

---

# **HVL Common Code Base Documentation**

*Release 0.10.0*

**Mikolaj Rybiński, David Graber, Henrik Menne, Alise Chachereau,**

**Jan 19, 2022**



## CONTENTS:

<b>1</b>	<b>HVL Common Code Base</b>	<b>1</b>
1.1	Features . . . . .	1
1.2	Documentation . . . . .	5
1.3	Credits . . . . .	5
<b>2</b>	<b>Installation</b>	<b>7</b>
2.1	Stable release . . . . .	7
2.2	From sources . . . . .	7
2.3	Additional system libraries . . . . .	8
<b>3</b>	<b>Usage</b>	<b>9</b>
<b>4</b>	<b>API Documentation</b>	<b>11</b>
4.1	hvl_ccb . . . . .	11
<b>5</b>	<b>Contributing</b>	<b>193</b>
5.1	Types of Contributions . . . . .	193
5.2	Get Started! . . . . .	194
5.3	Merge Request Guidelines . . . . .	195
5.4	Tips . . . . .	195
5.5	Deploying . . . . .	196
<b>6</b>	<b>Credits</b>	<b>197</b>
6.1	Maintainers . . . . .	197
6.2	Authors . . . . .	197
6.3	Contributors . . . . .	197
<b>7</b>	<b>History</b>	<b>199</b>
7.1	0.10.0 (2022-01-17) . . . . .	199
7.2	0.9.0 (2022-01-07) . . . . .	200
7.3	0.8.5 (2021-11-05) . . . . .	200
7.4	0.8.4 (2021-10-22) . . . . .	200
7.5	0.8.3 (2021-09-27) . . . . .	201
7.6	0.8.2 (2021-08-27) . . . . .	201
7.7	0.8.1 (2021-08-13) . . . . .	201
7.8	0.8.0 (2021-07-02) . . . . .	201
7.9	0.7.1 (2021-06-04) . . . . .	201
7.10	0.7.0 (2021-05-25) . . . . .	202
7.11	0.6.1 (2021-05-08) . . . . .	202
7.12	0.6.0 (2021-04-23) . . . . .	202
7.13	0.5.0 (2020-11-11) . . . . .	203

7.14	0.4.0 (2020-07-16)	204
7.15	0.3.5 (2020-02-18)	204
7.16	0.3.4 (2019-12-20)	204
7.17	0.3.3 (2019-05-08)	205
7.18	0.3.2 (2019-05-08)	205
7.19	0.3.1 (2019-05-02)	205
7.20	0.3 (2019-05-02)	205
7.21	0.2.1 (2019-04-01)	205
7.22	0.2.0 (2019-03-31)	205
7.23	0.1.0 (2019-02-06)	206
<b>8</b>	<b>Indices and tables</b>	<b>207</b>
	<b>Python Module Index</b>	<b>209</b>
	<b>Index</b>	<b>211</b>

## HVL COMMON CODE BASE

Python common code base (CCB) to control devices, which are used in high-voltage research. All implemented devices are used and tested in the High Voltage Laboratory (HVL) of the Federal Institute of Technology Zurich (ETH Zurich).

- Free software: GNU General Public License v3
- Copyright (c) 2019-2022 ETH Zurich, SIS ID and HVL D-ITET

### 1.1 Features

For managing multi-device experiments instantiate the `ExperimentManager` utility class.

#### 1.1.1 Devices

The device wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, various settings and options, as well as start/stop methods. Currently wrappers are available to control the following devices:

Function/Type	Devices
Data acquisition	LabJack (T4, T7, T7-PRO; requires <a href="#">LJM Library</a> ) Pico Technology PT-104 Platinum Resistance Data Logger (requires <a href="#">PicoSDK/libusbpt104</a> )
Digital Delay Generator	Highland T560
Digital IO	LabJack (T4, T7, T7-PRO; requires <a href="#">LJM Library</a> )
Experiment control	HVL Cube with and without Power Inverter
Gas Analyser	MBW 973-SF6 gas dew point mirror analyzer Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges SST Luminox oxygen sensor
I2C host	TiePie (HS5, WS5; requires <a href="#">LibTiePie SDK</a> )
Laser	CryLaS pulsed laser CryLaS laser attenuator
Oscilloscope	Rhode & Schwarz RTO 1024 TiePie (HS5, HS6, WS5; requires <a href="#">LibTiePie SDK</a> )
Power supply	Elektro-Automatik PSI9000 FuG Elektronik Heinzinger PNC Technix capacitor charger
Stepper motor drive	Newport SMC100PP Schneider Electric ILS2T
Temperature control	Lauda PRO RP 245 E circulation thermostat
Waveform generator	TiePie (HS5, WS5; requires <a href="#">LibTiePie SDK</a> )

Each device uses at least one standardised communication protocol wrapper.

### **1.1.2 Communication protocols**

In `hvl_ccb` by “communication protocol” we mean different levels of communication standards, from the low level actual communication protocols like serial communication to application level interfaces like VISA TCP standard. There are also devices in `hvl_ccb` that use a dummy communication protocol; this is because these devices are build on proprietary manufacturer libraries that communicate with the corresponding devices, as in the case of TiePie or LabJack devices.

The communication protocol wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, as well as open/close and read/write/query methods. Currently, wrappers for the following communication protocols are available:

Communication protocol	Devices using
Modbus TCP	Schneider Electric ILS2T stepper motor drive
OPC UA	HVL Cube with and without Power Inverter
Serial	<p>CryLaS pulsed laser and laser attenuator</p> <p>FuG Elektronik power supply (e.g. capacitor charger HCK) using the Probus V protocol</p> <p>Heinzinger PNC power supply using Heinzinger Digital Interface I/II</p> <p>SST Luminox oxygen sensor</p> <p>MBW 973-SF6 gas dew point mirror analyzer</p> <p>Newport SMC100PP single axis driver for 2-phase stepper motors</p> <p>Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges</p> <p>Technix capacitor charger</p>
TCP	Lauda PRO RP 245 E circulation thermostat
Telnet	Technix capacitor charger
VISA TCP	<p>Elektro-Automatik PSI9000 DC power supply</p> <p>Rhode &amp; Schwarz RTO 1024 oscilloscope</p>
<i>propriety</i>	<p>LabJack (T4, T7, T7-PRO) devices, which communicate via <a href="#">LJM Library</a></p> <p>Pico Technology PT-104 Platinum Resistance Data Logger, which communicate via <a href="#">PicoSDK/libusbpt104</a></p> <p>TiePie (HS5, HS6, WS5) oscilloscopes, generators and I2C hosts, which communicate via <a href="#">LibTiePie SDK</a></p>

### 1.1.3 Sensor and Unit Conversion Utility

The Sensor and Unit Conversion Utility is a submodule that allows on the one hand a unified implementation of hardware-sensors and on the other hand provides a unified way to convert units. The utility can be used with single numbers (`int`, `float`) as well as array-like structures containing single numbers (`np.array()`, `list`, `dict`, `tuple`).

Currently the following sensors are implemented:

- LEM LT 4000S
- LMT 70A

The following unit conversion classes are implemented:

- Temperature (Kelvin, Celsius, Fahrenheit)
- Pressure (Pascal, Bar, Atmosphere, Psi, Torr, Millimeter Mercury)

## 1.2 Documentation

Note: if you're planning to contribute to the `hvl_ccb` project read the **Contributing** section in the HVL CCB documentation.

Do either:

- read [HVL CCB documentation at RTD](#),

or

- build and read HVL CCB documentation locally; install first [Graphviz](#) (make sure to have the `dot` command in the executable search path) and the Python build requirements for documentation:

```
$ pip install docs/requirements.txt
```

and then either on Windows in Git BASH run:

```
$ ./make.sh docs
```

or from any other shell with GNU Make installed run:

```
$ make docs
```

The target index HTML ("`docs/_build/html/index.html`") should open automatically in your Web browser.

## 1.3 Credits

This package was created with [Cookiecutter](#) and the [audreyr/cookiecutter-pypackage](#) project template.



## INSTALLATION

### 2.1 Stable release

To install HVL Common Code Base, run this command in your terminal:

```
$ pip install hvl_ccb
```

To install HVL Common Code Base with optional Python libraries that require manual installations of additional system libraries, you need to specify on installation extra requirements corresponding to these controllers. For instance, to install Python requirements for LabJack and TiePie devices, run:

```
$ pip install "hvl_ccb[tiapie,labjack]"
```

See below for the info about additional system libraries and the corresponding extra requirements.

To install all extra requirements run:

```
$ pip install "hvl_ccb[all]"
```

This is the preferred method to install HVL Common Code Base, as it will always install the most recent stable release. If you don't have `pip` installed, this [Python installation guide](#) can guide you through the process.

### 2.2 From sources

The sources for HVL Common Code Base can be downloaded from the [GitLab repo](#).

You can either clone the repository:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

Or download the `tarball`:

```
$ curl -OL https://gitlab.com/ethz_hvl/hvl_ccb/-/archive/master/hvl_ccb.tar.gz
```

Once you have a copy of the source, you can install it with:

```
$ pip install .
```

## 2.3 Additional system libraries

If you have installed *hvl\_ccb* with any of the extra features corresponding to device controllers, you must additionally install respective system library; these are:

Extra feature	Additional system library
labjack	<a href="#">LJM Library</a>
picotech	<a href="#">PicoSDK (Windows)</a> / <a href="#">libusbpt104 (Ubuntu/Debian)</a>
tiepie	<a href="#">LibTiePie SDK</a>

For more details on installation of the libraries see docstrings of the corresponding *hvl\_ccb* modules.

---

CHAPTER  
**THREE**

---

**USAGE**

To use HVL Common Code Base in a project:

```
import hvl_ccb
```



## API DOCUMENTATION


### 4.1 hvl\_ccb

#### 4.1.1 Subpackages

`hvl_ccb.comm`

Submodules

`hvl_ccb.comm.base`

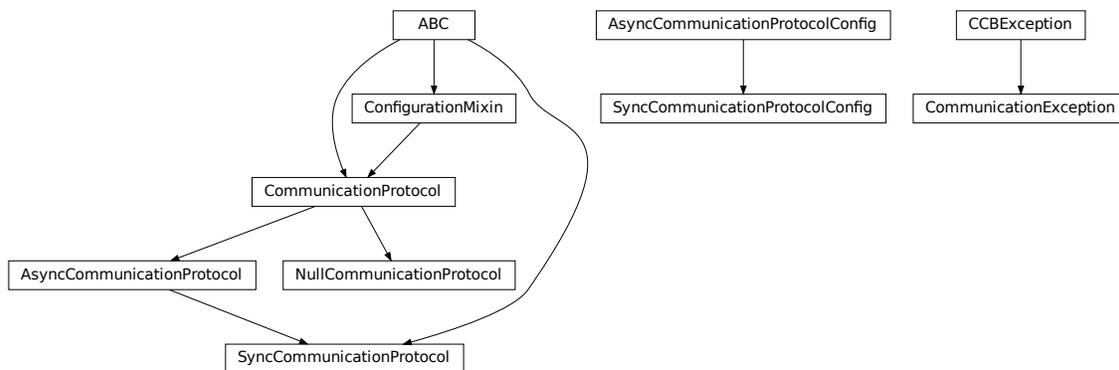

Module with base classes for communication protocols.

**class** `AsyncCommunicationProtocol`(*config*)

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Abstract base class for asynchronous communication protocols

**static** `config_cls()` → `Type[hvl_ccb.comm.base.AsyncCommunicationProtocolConfig]`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**read()** → `str`

Read a single line of text as *str* from the communication.

**Returns** text as *str* including the terminator, which can also be empty ""

**read\_all**(*n\_attempts\_max*: *Optional[int] = None*, *attempt\_interval\_sec*: *Optional[Union[int, float]] = None*) → *Optional[str]*

Read all lines of text from the connection till nothing is left to read.

**Parameters**

- **n\_attempts\_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt\_interval\_sec** – time between the reading attempts

**Returns** A multi-line *str* including the terminator internally

**abstract read\_bytes**() → *bytes*

Read a single line as *bytes* from the communication.

This method uses *self.access\_lock* to ensure thread-safety.

**Returns** a single line as *bytes* containing the terminator, which can also be empty b''''

**read\_nonempty**(*n\_attempts\_max*: *Optional[int] = None*, *attempt\_interval\_sec*: *Optional[Union[int, float]] = None*) → *Optional[str]*

Try to read a non-empty single line of text as *str* from the communication. If the host does not reply or reply with white space only, it will return *None*.

**Returns** a non-empty text as a *str* or *None* in case of an empty string

**Parameters**

- **n\_attempts\_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt\_interval\_sec** – time between the reading attempts

**read\_text**() → *str*

Read one line of text from the serial port. The input buffer may hold additional data afterwards, since only one line is read.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

**Returns** String read from the serial port; '' if there was nothing to read.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**read\_text\_nonempty**(*n\_attempts\_max*: *Optional[int] = None*, *attempt\_interval\_sec*: *Optional[Union[int, float]] = None*) → *Optional[str]*

Reads from the serial port, until a non-empty line is found, or the number of attempts is exceeded.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Attention: in contrast to *read\_text*, the returned answer will be stripped of a whitespace newline terminator at the end, if such terminator is set in the initial configuration (default).

**Parameters**

- **n\_attempts\_max** – maximum number of read attempts
- **attempt\_interval\_sec** – time between the reading attempts

**Returns** String read from the serial port; '' if number of attempts is exceeded or serial port is not opened.

**write**(*text*: *str*)

Write text as *str* to the communication.

**Parameters** **text** – text as a *str* to be written

**abstract write\_bytes**(*data: bytes*) → int  
Write data as *bytes* to the communication.

This method uses *self.access\_lock* to ensure thread-safety.

**Parameters** **data** – data as *bytes*-string to be written

**Returns** number of bytes written

**write\_text**(*text: str*)

Write text to the serial port. The text is encoded and terminated by the configured terminator.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

**Parameters** **text** – Text to send to the port.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

```
class AsyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                     encoding_error_handling: str = 'strict',
                                     wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                     default_n_attempts_read_text_nonempty: int = 10)
```

Bases: object

Base configuration data class for asynchronous communication protocols

**clean\_values**()

**default\_n\_attempts\_read\_text\_nonempty: int = 10**  
default number of attempts to read a non-empty text

**encoding: str = 'utf-8'**  
Standard encoding of the connection. Typically this is *utf-8*, but can also be *latin-1* or something from here: <https://docs.python.org/3/library/codecs.html#standard-encodings>

**encoding\_error\_handling: str = 'strict'**  
Encoding error handling scheme as defined here: <https://docs.python.org/3/library/codecs.html#error-handlers> By default strict error handling that raises *UnicodeError*.

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys**() → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod** `required_keys()` → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**terminator:** bytes = b'\r\n'

The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations. This defines the end of a single line.

**wait\_sec\_read\_text\_nonempty:** Union[int, float] = 0.5

time to wait between attempts of reading a non-empty text

**exception** `CommunicationException`

Bases: `hvl_ccb.exception.CCBException`

**class** `CommunicationProtocol(config)`

Bases: `hvl_ccb.configuration.ConfigurationMixin`, `abc.ABC`

Communication protocol abstract base class.

Specifies the methods to implement for communication protocol, as well as implements some default settings and checks.

**access\_lock**

Access lock to use with context manager when accessing the communication protocol (thread safety)

**abstract** `close()`

Close the communication protocol

**abstract** `open()`

Open communication protocol

**class** `NullCommunicationProtocol(config)`

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Communication protocol that does nothing.

**close()** → None

Void close function.

**static** `config_cls()` → Type[`hvl_ccb.configuration.EmptyConfig`]

Empty configuration

**Returns** EmptyConfig

**open()** → None

Void open function.

**class** `SyncCommunicationProtocol(config)`

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocol`, `abc.ABC`

Abstract base class for synchronous communication protocols with `query()`

**static** `config_cls()` → Type[`hvl_ccb.comm.base.SyncCommunicationProtocolConfig`]

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query**(*command: str*) → Optional[str]

Send a command to the interface and handle the status message. Eventually raises an exception.


**Parameters** `command` – Command to send

**Returns** Answer from the interface, which can be None instead of an empty reply

```
class SyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                     encoding_error_handling: str = 'strict',
                                     wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                     default_n_attempts_read_text_nonempty: int = 10)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

## hvl\_ccb.comm.labjack\_ljm

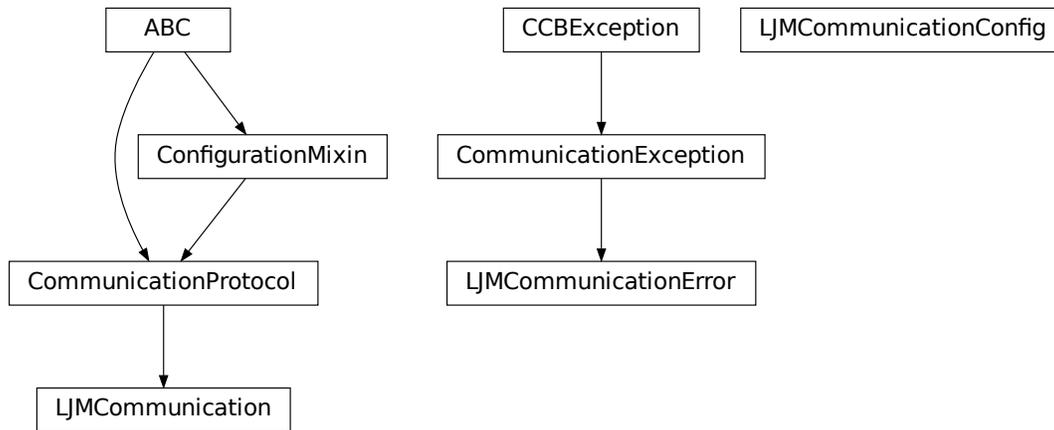

Communication protocol for LabJack using the LJM Library. Originally developed and tested for LabJack T7-PRO.

Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

```
class LJMCommunication(configuration)
```

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Communication protocol implementing the LabJack LJM Library Python wrapper.

**close()** → *None*

Close the communication port.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property is\_open: bool**

Flag indicating if the communication port is open.

**Returns** *True* if the port is open, otherwise *False*

**open()** → *None*

Open the communication port.

**read\_name(\*names: str, return\_num\_type: typing.Type[numbers.Real] = <class 'float'>) → Union[numbers.Real, Sequence[numbers.Real]]**

Read one or more input numeric values by name.

**Parameters**

- **names** – one or more names to read out from the LabJack
- **return\_num\_type** – optional numeric type specification for return values; by default *float*.

**Returns** answer of the LabJack, either single number or multiple numbers in a sequence, respectively, when one or multiple names to read were given

**Raises** **TypeError** – if read value of type not compatible with *return\_num\_type*

**write\_name**(*name: str, value: numbers.Real*) → *None*

Write one value to a named output.

**Parameters**

- **name** – String or with name of LabJack IO
- **value** – is the value to write to the named IO port

**write\_names**(*name\_value\_dict: Dict[str, numbers.Real]*) → *None*

Write more than one value at once to named outputs.

**Parameters** **name\_value\_dict** – is a dictionary with string names of LabJack IO as keys and corresponding numeric values

```
class LJMCommunicationConfig(device_type: Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY',  
                             connection_type: Union[str,  
                             hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType] =  
                             'ANY', identifier: str = 'ANY')
```

Bases: object

Configuration dataclass for *LJMCommunication*.

```
class ConnectionType(value=<no_arg>, names=None, module=None, type=None, start=1,  
                    boundary=None)
```

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

LabJack connection type.

**ANY** = 1

**ETHERNET** = 4

**TCP** = 3

**USB** = 2

**WIFI** = 5

```
class DeviceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

LabJack device types.

Can be also looked up by ambiguous Product ID (*p\_id*) or by instance name: ``python LabJackDeviceType(4) is LabJackDeviceType('T4')``

**ANY** = 1

**T4** = 2

**T7** = 3

**T7\_PRO** = 4

```
classmethod get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,  
List[hvl_ccb._dev.labjack.DeviceType]]
```

Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

**Parameters** **p\_id** – Product ID of a LabJack device

**Returns** Instance or list of instances of *LabJackDeviceType*

**Raises** **ValueError** – when Product ID is unknown

```
clean_values() → None
```

Performs value checks on *device\_type* and *connection\_type*.

```
connection_type: Union[str,  
hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType] = 'ANY'
```

Can be either string or of enum *ConnectionType*.

```
device_type: Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY'
```

Can be either string 'ANY', 'T7\_PRO', 'T7', 'T4', or of enum *DeviceType*.

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
identifier: str = 'ANY'
```

The identifier specifies information for the connection to be used. This can be an IP address, serial number, or device name. See the LabJack docs ( <https://labjack.com/support/software/api/ljm/function-reference/ljmopens/identifier-parameter>) for more information.

```
is_configdataclass = True
```

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

```
classmethod optional_defaults() → Dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


**Returns** a list of strings containing all required keys.

```
exception LJMCommunicationError
```

Bases: *hvl\_ccb.comm.base.CommunicationException*

Errors coming from LJMCommunication.

## hvl\_ccb.comm.modbus\_tcp

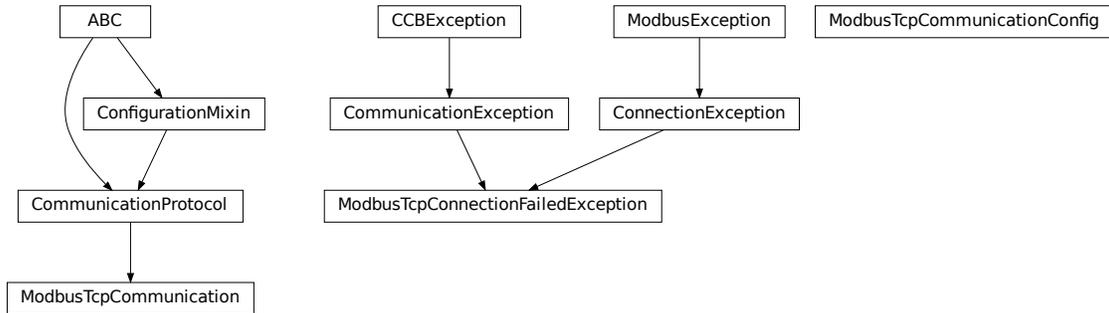

Communication protocol for modbus TCP ports. Makes use of the [pymodbus](#) library.

**class** `ModbusTcpCommunication(configuration)`

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Implements the Communication Protocol for modbus TCP.

**close()**

Close the Modbus TCP connection.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**open()** → *None*

Open the Modbus TCP connection.

**Raises** `ModbusTcpConnectionFailedException` – if the connection fails.

**read\_holding\_registers(address: int, count: int)** → List[int]

Read specified number of register starting with given address and return the values from each register.

**Parameters**

- **address** – address of the first register
- **count** – count of registers to read

**Returns** list of *int* values

**read\_input\_registers(address: int, count: int)** → List[int]

Read specified number of register starting with given address and return the values from each register in a list.

**Parameters**

- **address** – address of the first register
- **count** – count of registers to read

**Returns** list of *int* values

**write\_registers(address: int, values: Union[List[int], int])**

Write values from the specified address forward.

**Parameters**

- **address** – address of the first register
- **values** – list with all values

**class ModbusTcpCommunicationConfig**(*host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], unit: int, port: int = 502*)

Bases: object

Configuration dataclass for *ModbusTcpCommunication*.

**clean\_values()**

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**host:** Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]

**is\_configdataclass** = True

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port:** int = 502

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


**unit:** int

**exception ModbusTcpConnectionFailedException**(*string=""*)

Bases: pymodbus.exceptions.ConnectionException, *hvl\_ccb.comm.base.CommunicationException*

Exception raised when the connection failed.

## hvl\_ccb.comm.opc

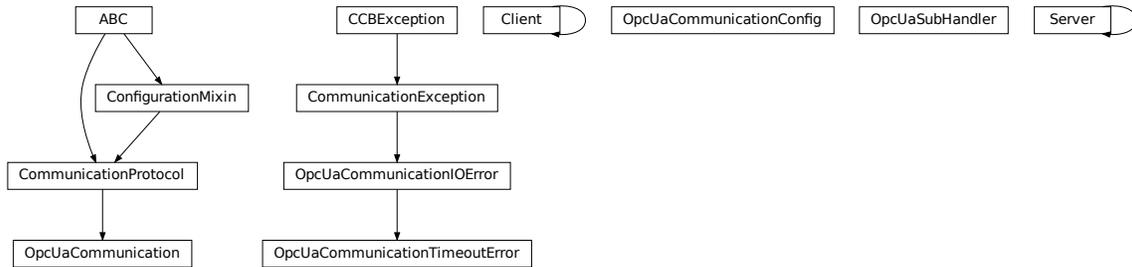

Communication protocol implementing an OPC UA connection. This protocol is used to interface with the “Super-cube” PLC from Siemens.

```
class Client(url: str, timeout: int = 4)
```

Bases: `asyncua.sync.Client`

```
disconnect()
```

```
get_objects_node()
```

Get Objects node of client. Returns a Node object.

```
property is_open
```

```
send_hello(*args, **kwargs)
```

```
class OpcUaCommunication(config)
```

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Communication protocol implementing an OPC UA connection. Makes use of the package python-opcua.

```
close() → None
```

Close the connection to the OPC UA server.

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
init_monitored_nodes(node_id: Union[object, Iterable], ns_index: int) → None
```

Initialize monitored nodes.

#### Parameters

- **node\_id** – one or more strings of node IDs; node IDs are always casted via `str()` method here, hence do not have to be strictly string objects.
- **ns\_index** – the namespace index the nodes belong to.

**Raises** `OpcUaCommunicationIOError` – when protocol was not opened or can’t communicate with a OPC UA server

```
property is_open: bool
```

Flag indicating if the communication port is open.

**Returns** `True` if the port is open, otherwise `False`

```
open() → None
```

Open the communication to the OPC UA server.

Raises *OpcUaCommunicationIOError* – when communication port cannot be opened.

**read**(*node\_id, ns\_index*)

Read a value from a node with id and namespace index.

**Parameters**

- **node\_id** – the ID of the node to read the value from
- **ns\_index** – the namespace index of the node

**Returns** the value of the node object.

Raises *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

**write**(*node\_id, ns\_index, value*) → *None*

Write a value to a node with name *name*.

**Parameters**

- **node\_id** – the id of the node to write the value to.
- **ns\_index** – the namespace index of the node.
- **value** – the value to write.

Raises *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

```
class OpcUaCommunicationConfig(host: typing.Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],
                               endpoint_name: str, port: int = 4840, sub_handler:
                               hvl_ccb.comm.opc.OpcUaSubHandler =
                               <hvl_ccb.comm.opc.OpcUaSubHandler object>, update_parameter:
                               asyncua.ua.uaprotoocol_auto.CreateSubscriptionParameters =
                               CreateSubscriptionParameters(RequestedPublishingInterval=1000,
                                                             RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22,
                                                             MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0),
                               wait_timeout_retry_sec: typing.Union[int, float] = 1,
                               max_timeout_retry_nr: int = 5)
```

Bases: object

Configuration dataclass for OPC UA Communciation.

**clean\_values**()

**endpoint\_name: str**

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]**

Hostname or IP-Address of the OPC UA server.

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**max\_timeout\_retry\_nr: int = 5**

Maximal number of call re-tries on underlying OPC UA client timeout error

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 4840**

Port of the OPC UA server to connect to.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**sub\_handler: hv1\_ccb.comm.opc.OpcUaSubHandler = <hv1\_ccb.comm.opc.OpcUaSubHandler object>**

object to use for handling subscriptions.

**update\_parameter: asyncua.ua.uaprotocol\_auto.CreateSubscriptionParameters = CreateSubscriptionParameters(RequestedPublishingInterval=1000, RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22, MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0)**

Values are given as a *ua.CreateSubscriptionParameters* as these parameters are requested by the OPC server. Other values will lead to an automatic revision of the parameters and a warning in the opc-logger, cf. MR !173

**wait\_timeout\_retry\_sec: Union[int, float] = 1**

Wait time between re-trying calls on underlying OPC UA client timeout error

**exception OpcUaCommunicationIOError**

Bases: OSError, *hv1\_ccb.comm.base.CommunicationException*

OPC-UA communication I/O error.

**exception OpcUaCommunicationTimeoutError**

Bases: *hv1\_ccb.comm.opc.OpcUaCommunicationIOError*

OPC-UA communication timeout error.

**class OpcUaSubHandler**

Bases: object

Base class for subscription handling of OPC events and data change events. Override methods from this class to add own handling capabilities.

To receive events from server for a subscription data\_change and event methods are called directly from receiving thread. Do not do expensive, slow or network operation there. Create another thread if you need to do such a thing.

**datachange\_notification**(node, val, data)


**event\_notification**(event)

**class Server**(shelf\_file=None, tloop=None)

Bases: *asyncua.sync.Server*

**get\_objects\_node()**

Get Objects node of server. Returns a Node object.

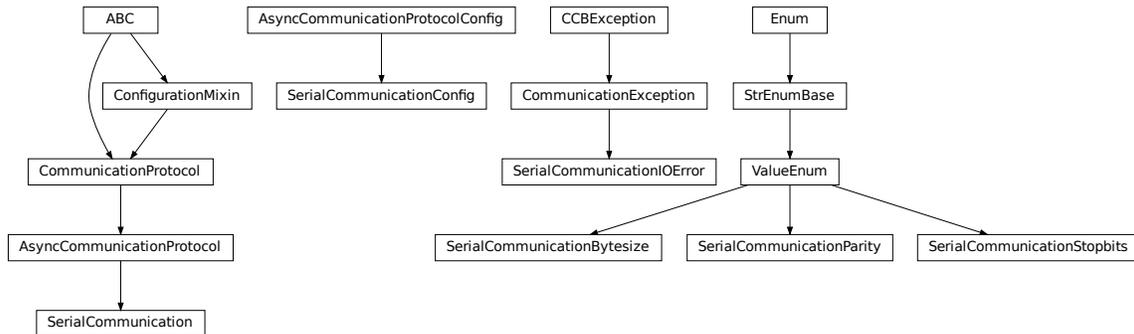**hvl\_ccb.comm.serial**

Communication protocol for serial ports. Makes use of the `pySerial` library.

**class SerialCommunication(configuration)**

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocol`

Implements the Communication Protocol for serial ports.

**close()**

Close the serial connection.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property is\_open: bool**

Flag indicating if the serial port is open.

**Returns** `True` if the serial port is open, otherwise `False`

**open()**

Open the serial connection.

**Raises** `SerialCommunicationIOError` – when communication port cannot be opened.

**read\_bytes() → bytes**

Read the bytes from the serial port till the terminator is found. The input buffer may hold additional lines afterwards.

This method uses `self.access_lock` to ensure thread-safety.

**Returns** Bytes read from the serial port; `b''` if there was nothing to read.

**Raises** `SerialCommunicationIOError` – when communication port is not opened

**read\_single\_bytes(size: int = 1) → bytes**

Read the specified number of bytes from the serial port. The input buffer may hold additional data afterwards.

**Returns** Bytes read from the serial port; `b''` if there was nothing to read.

**write\_bytes**(*data: bytes*) → int  
Write bytes to the serial port.

This method uses *self.access\_lock* to ensure thread-safety.

**Parameters** *data* – data to write to the serial port

**Returns** number of bytes written

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**class** *SerialCommunicationBytesize*(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*,  
*boundary=None*)

Bases: *hvl\_ccb.utils.enum.ValueEnum*

Serial communication bytesize.

**EIGHTBITS** = 8

**FIVEBITS** = 5

**SEVENBITS** = 7

**SIXBITS** = 6

**class** *SerialCommunicationConfig*(*terminator: bytes = b'\r\n'*, *encoding: str = 'utf-8'*,  
*encoding\_error\_handling: str = 'strict'*, *wait\_sec\_read\_text\_nonempty:*  
*Union[int, float] = 0.5*, *default\_n\_attempts\_read\_text\_nonempty: int = 10*,  
*port: Optional[str] = None*, *baudrate: int = 9600*, *parity: Union[str,*  
*hvl\_ccb.comm.serial.SerialCommunicationParity] =*  
*SerialCommunicationParity.NONE*, *stopbits: Union[int, float,*  
*hvl\_ccb.comm.serial.SerialCommunicationStopbits] =*  
*SerialCommunicationStopbits.ONE*, *bytesize: Union[int,*  
*hvl\_ccb.comm.serial.SerialCommunicationBytesize] =*  
*SerialCommunicationBytesize.EIGHTBITS*, *timeout: Union[int, float] = 2*)

Bases: *hvl\_ccb.comm.base.AsyncCommunicationProtocolConfig*

Configuration dataclass for *SerialCommunication*.

**Bytesize**

alias of *hvl\_ccb.comm.serial.SerialCommunicationBytesize*

**Parity**

alias of *hvl\_ccb.comm.serial.SerialCommunicationParity*

**Stopbits**

alias of *hvl\_ccb.comm.serial.SerialCommunicationStopbits*

**baudrate: int = 9600**

Baudrate of the serial port

**bytesize: Union[int, *hvl\_ccb.comm.serial.SerialCommunicationBytesize*] = 8**

Size of a byte, 5 to 8

**clean\_values()**

**create\_serial\_port()** → *serial.serialposix.Serial*

Create a serial port instance according to specification in this configuration

**Returns** Closed serial port instance

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** Union[str, *hvl\_ccb.comm.serial.SerialCommunicationParity*] = 'N'

Parity to be used for the connection.

**port:** Optional[str] = None

Port is a string referring to a COM-port (e.g. 'COM3') or a URL. The full list of capabilities is found on [the pyserial documentation](#).

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** Union[int, float, *hvl\_ccb.comm.serial.SerialCommunicationStopbits*] = 1

Stopbits setting, can be 1, 1.5 or 2.

**terminator\_str()** → str

**timeout:** Union[int, float] = 2

Timeout in seconds for the serial port

**exception SerialCommunicationIOError**

Bases: OSError, *hvl\_ccb.comm.base.CommunicationException*

Serial communication related I/O errors.

**class SerialCommunicationParity**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl\_ccb.utils.enum.ValueEnum*

Serial communication parity.

**EVEN** = 'E'

**MARK** = 'M'

**NAMES** = {'E': 'Even', 'M': 'Mark', 'N': 'None', 'O': 'Odd', 'S': 'Space'}

**NONE** = 'N'

**ODD** = 'O'

**SPACE** = 'S'

```
class SerialCommunicationStopbits(value=<no_arg>, names=None, module=None, type=None, start=1,
                                boundary=None)
```

Bases: *hvl\_ccb.utils.enum.ValueEnum*


Serial communication stopbits.

**ONE** = 1

**ONE\_POINT\_FIVE** = 1.5

**TWO** = 2

### hvl\_ccb.comm.tcp

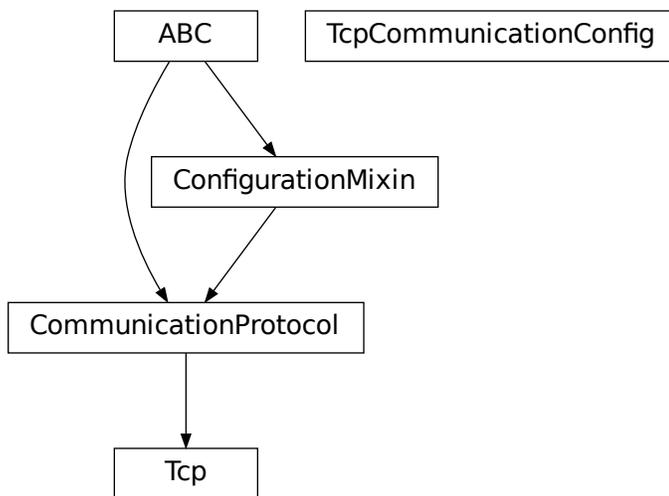

TCP communication protocol.

Makes use of the socket library.

```
class Tcp(configuration)
```

Bases: *hvl\_ccb.comm.base.CommunicationProtocol*

Tcp Communication Protocol.

**close()** → *None*

Close TCP connection.

**static config\_cls()** → *Type[hvl\_ccb.comm.tcp.TcpCommunicationConfig]*

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**open()** → *None*

Open TCP connection.

**read()** → str

TCP read function :return: information read from TCP buffer formatted as string

**write**(*command: str = ""*) → *None*

TCP write function :param *command*: command string to be sent :return: none

**class TcpCommunicationConfig**(*host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]*, *port: int = 54321*, *bufsize: int = 1024*)

Bases: object

Configuration dataclass for TcpCommunication.

**bufsize: int = 1024**

**clean\_values()**

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]**

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.


**port: int = 54321**

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

## hvl\_ccb.comm.telnet

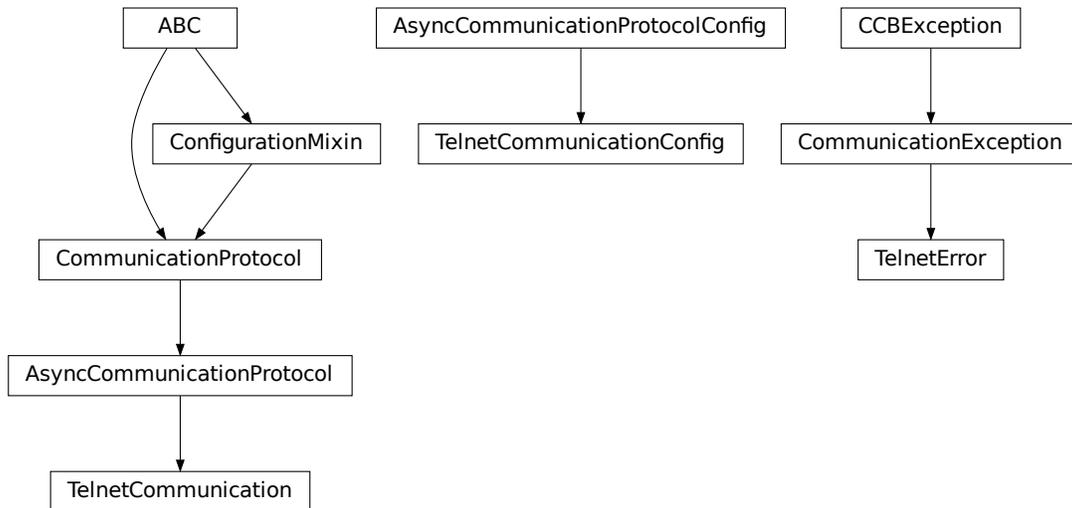

Communication protocol for telnet. Makes use of the [telnetlib](#) library.

**class TelnetCommunication**(*configuration*)

Bases: [hvl\\_ccb.comm.base.AsyncCommunicationProtocol](#)

Implements the Communication Protocol for telnet.

**close()**

Close the telnet connection unless it is not closed.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property is\_open: bool**

Is the connection open?

**Returns** True for an open connection

**open()**

Open the telnet connection unless it is not yet opened.

**read\_bytes()** → bytes

Read data as *bytes* from the telnet connection.

**Returns** data from telnet connection

**Raises** [TelnetError](#) – when connection is not open, raises an Error during the communication

**write\_bytes(data: bytes)**

Write the data as *bytes* to the telnet connection.

**Parameters data** – Data to be sent.

**Raises** [TelnetError](#) – when connection is not open, raises an Error during the communication

```
class TelnetCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,
                                host: Optional[Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]]
                                = None, port: int = 0, timeout: Union[int, float] = 0.2)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

Configuration dataclass for `TelnetCommunication`.

**clean\_values()**

**create\_telnet()** → `Optional[telnetlib.Telnet]`

Create a telnet client :return: Opened Telnet object or None if connection is not possible

**force\_value(*fieldname, value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**host: Optional[Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]] = None**

Host to connect to can be localhost or

**classmethod keys()** → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → `Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 0**

Port at which the host is listening

**classmethod required\_keys()** → `Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**timeout: Union[int, float] = 0.2**


Timeout for reading a line

**exception TelnetError**

Bases: `OSError`, `hvl_ccb.comm.base.CommunicationException`

Telnet communication related errors.

## hvl\_ccb.comm.visa

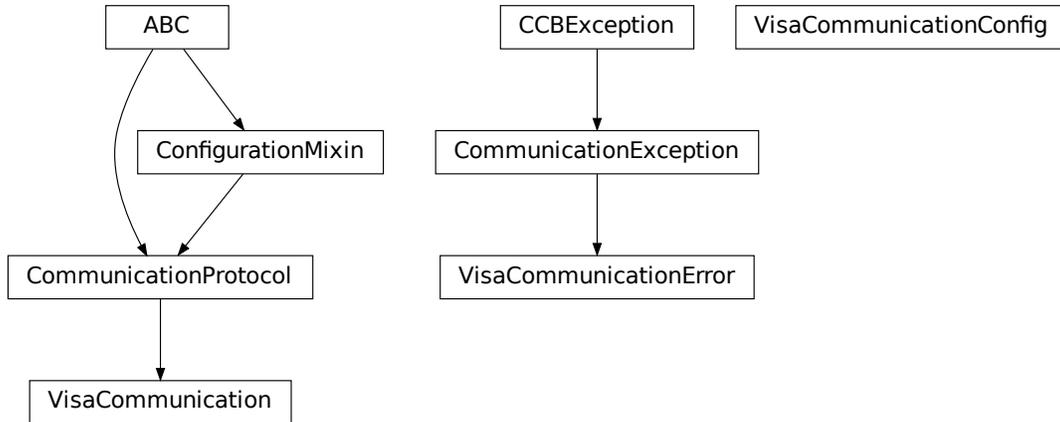

Communication protocol for VISA. Makes use of the pyvisa library. The backend can be NI-Visa or pyvisa-py.

Information on how to install a VISA backend can be found here: [https://pyvisa.readthedocs.io/en/master/getting\\_nivisa.html](https://pyvisa.readthedocs.io/en/master/getting_nivisa.html)

So far only TCPIP SOCKET and TCPIP INSTR interfaces are supported.

**class VisaCommunication**(*configuration*)

Bases: *hvl\_ccb.comm.base.CommunicationProtocol*

Implements the Communication Protocol for VISA / SCPI.

**MULTI\_COMMANDS\_MAX** = 5

The maximum of commands that can be sent in one round is 5 according to the VISA standard.

**MULTI\_COMMANDS\_SEPARATOR** = ';'

The character to separate two commands is ; according to the VISA standard.

**WAIT\_AFTER\_WRITE** = 0.08

Small pause in seconds to wait after write operations, allowing devices to really do what we tell them before continuing with further tasks.

**close()** → *None*

Close the VISA connection and invalidates the handle.

**static config\_cls()** → *Type[hvl\_ccb.comm.visa.VisaCommunicationConfig]*

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**open()** → *None*

Open the VISA connection and create the resource.

**query**(\**commands: str*) → *Union[str, Tuple[str, ...]]*

A combination of write(message) and read.

**Parameters** **commands** – list of commands

**Returns** list of values

**Raises** *VisaCommunicationError* – when connection was not started, or when trying to issue too many commands at once.

**spoll()** → int

Execute serial poll on the device. Reads the status byte register STB. This is a fast function that can be executed periodically in a polling fashion.

**Returns** integer representation of the status byte

**Raises** *VisaCommunicationError* – when connection was not started

**write(\*commands: str)** → None

Write commands. No answer is read or expected.

**Parameters** **commands** – one or more commands to send

**Raises** *VisaCommunicationError* – when connection was not started

```
class VisaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],
                              interface_type: Union[str,
                                                    hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType], board: int =
0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800,
                              open_timeout: int = 1000, write_termination: str = '\n', read_termination: str
                              = '\n', visa_backend: str = "")
```

Bases: object

*VisaCommunication* configuration dataclass.

```
class InterfaceType(value=<no_arg>, names=None, module=None, type=None, start=1,
                    boundary=None)
```

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

Supported VISA Interface types.

**TCPIP\_INSTR** = 2

VXI-11 protocol

**TCPIP\_SOCKET** = 1

VISA-RAW protocol

**address**(host: str, port: Optional[int] = None, board: Optional[int] = None) → str

Address string specific to the VISA interface type.

**Parameters**

- **host** – host IP address
- **port** – optional TCP port
- **board** – optional board number

**Returns** address string

**property address:** str

Address string depending on the VISA protocol's configuration.

**Returns** address string corresponding to current configuration

**board:** int = 0

Board number is typically 0 and comes from old bus systems.

**chunk\_size:** int = 204800

Chunk size is the allocated memory for read operations. The standard is 20kB, and is increased per default here to 200kB. It is specified in bytes.

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host**: Union[str, `ipaddress.IPv4Address`, `ipaddress.IPv6Address`]

**interface\_type**: Union[str, `hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType`]

Interface type of the VISA connection, being one of *InterfaceType*.

**is\_configdataclass** = True

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**open\_timeout**: int = 1000

Timeout for opening the connection, in milli seconds.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port**: int = 5025

TCP port, standard is 5025.

**read\_termination**: str = '\n'

Read termination character.

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**timeout**: int = 5000

Timeout for commands in milli seconds.

**visa\_backend**: str = ''

Specifies the path to the library to be used with PyVISA as a backend. Defaults to None, which is NI-VISA (if installed), or pyvisa-py (if NI-VISA is not found). To force the use of pyvisa-py, specify '@py' here.

**write\_termination**: str = '\n'

Write termination character.

**exception VisaCommunicationError**

Bases: OSError, `hvl_ccb.comm.base.CommunicationException`

Base class for VisaCommunication errors.

## Module contents

Communication protocols subpackage.


`hvl_ccb.dev`

## Subpackages

`hvl_ccb.dev.cube`

## Submodules

`hvl_ccb.dev.cube.base`

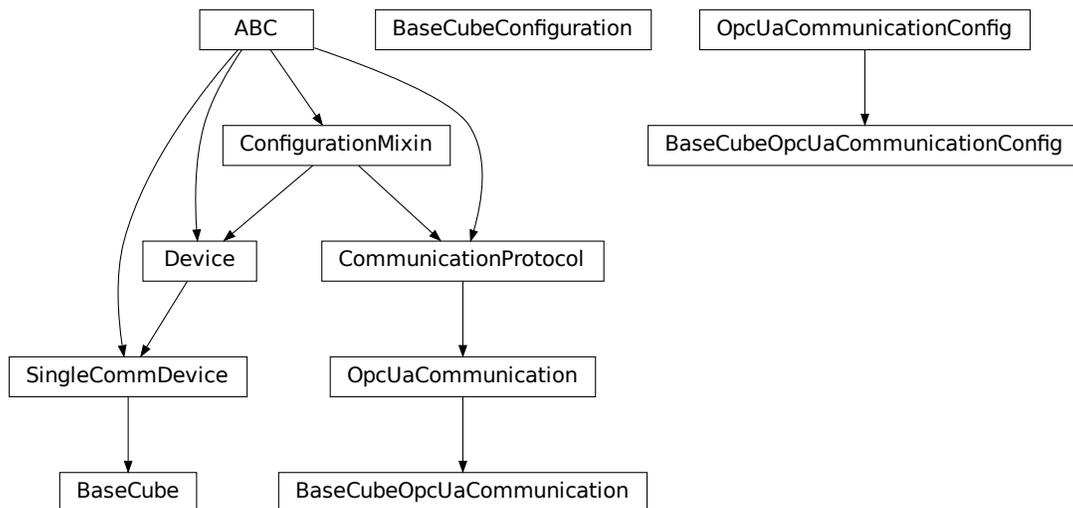

Classes for the BaseCube device.

```
class BaseCube(com, dev_config=None)
```

Bases: `hvl_ccb.dev.base.SingleCommDevice`

Base class for Cube variants.

**OPC\_MAX\_YEAR** = 2089

**OPC\_MIN\_YEAR** = 1990

**active\_alarms**(*human\_readable*: *bool* = True) → List[Union[int, str]]

Displays all active alarms / messages.

**Parameters** *human\_readable* – True for human readable message, False for corresponding integer

**Returns** list with active alarms

**property breakdown\_detection\_active: bool**

Get the state of the breakdown detection functionality. Returns True if it is enabled, False otherwise.

**Returns** state of the breakdown detection functionality

**breakdown\_detection\_reset()** → *None*

Reset the breakdown detection circuitry so that it is ready to detect breakdowns again.

**property breakdown\_detection\_triggered: bool**

See if breakdown detection unit has been triggered. Returns True if it is triggered, False otherwise.

**Returns** trigger status of the breakdown detection unit

**property cee16\_socket**

Read the on-state of the IEC CEE16 three-phase power socket.

**Returns** the on-state of the CEE16 power socket

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**classmethod datetime\_to\_opc(time\_dt: datetime.datetime)** → List[int]

Converts python datetime format into opc format (list of 8 integers) as defined in the following link: <https://support.industry.siemens.com/cs/mdm/109798671?c=133950752267&lc=de-WW> Each byte corresponds to one list entry. [yy, MM, dd, hh, mm, ss, milliseconds, weekday] Milliseconds and Weekday are not used, as this precision / information is not needed. The conversion of the numbers is special. Each decimal number is treated as it would be a hex-number and then converted back to decimal. This is tested with the used PLC in the BaseCube. yy: 0 to 99 (0 -> 2000, 89 -> 2089, 90 -> 1990, 99 -> 1999) MM: 1 to 12 dd: 1 to 31 hh: 0 to 23 mm: 0 to 59 ss: 0 to 59

**Parameters time\_dt** – time to be converted

**Returns** time in opc list format

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**display\_message\_board()** → *None*

Display 15 newest messages

**display\_status\_board()** → *None*

Display status board.

**door\_1\_status**

Get the status of a safety fence door. See constants.DoorStatus for possible returned door statuses.

**door\_2\_status**

Get the status of a safety fence door. See constants.DoorStatus for possible returned door statuses.

**door\_3\_status**

Get the status of a safety fence door. See constants.DoorStatus for possible returned door statuses.

**earthing\_rod\_1\_status**

Get the status of a earthing rod. See constants.EarthingRodStatus for possible returned earthing rod statuses.

**earthing\_rod\_2\_status**

Get the status of a earthing rod. See constants.EarthingRodStatus for possible returned earthing rod statuses.

**earthing\_rod\_3\_status**

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

**property operate: Optional[bool]**

Indicates if 'operate' is activated. 'operate' means locket safety circuit, red lamps, high voltage on and locked safety switches.

**Returns** *True* if operate is activated (`RedOperate`), *False* if ready is deactivated (`RedReady`), *None* otherwise

**quit\_error()** → *None*

Quits errors that are active on the Cube.

**read(node\_id: str)**

Local wrapper for the OPC UA communication protocol read method.

**Parameters** `node_id` – the id of the node to read.

**Returns** the value of the variable

**property ready: Optional[bool]**

Indicates if 'ready' is activated. 'ready' means locket safety circuit, red lamps, but high voltage still off.

**Returns** *True* if ready is activated (`RedReady`), *False* if ready is deactivated (`GreenReady`), *None* otherwise

**set\_message\_board(msgs: List[str], display\_board: bool = True) → None**

Fills messages into message board that display that 15 newest messages with a timestamp.

**Parameters**

- `msgs` – list of strings
- `display_board` – display 15 newest messages if *True* (default)

**Raises** `ValueError` – if there are too many messages or the positions indices are invalid.

**set\_status\_board(msgs: List[str], pos: Optional[List[int]] = None, clear\_board: bool = True, display\_board: bool = True) → None**

Sets and displays a status board. The messages and the position of the message can be defined.

**Parameters**

- `msgs` – list of strings
- `pos` – list of integers [0...14]
- `clear_board` – clear unspecified lines if *True* (default), keep otherwise
- `display_board` – display new status board if *True* (default)

**Raises** `ValueError` – if there are too many messages or the positions indices are invalid.

**start()** → *None*

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

**property status: hvl\_ccb.dev.cube.constants.SafetyStatus**

Get the safety circuit status of the Cube. This methods is for the user.

**Returns** the safety status of the Cube's state machine.

**stop()** → *None*

Stop the Cube device. Deactivates the remote control and closes the communication protocol.

Raises *CubeStopError* – when the cube is not in the correct status to stop the operation

**t13\_socket\_1**

Set and get the state of a SEV T13 power socket.

**t13\_socket\_2**

Set and get the state of a SEV T13 power socket.

**t13\_socket\_3**

Set and get the state of a SEV T13 power socket.

**write**(*node\_id*, *value*) → *None*

Local wrapper for the OPC UA communication protocol write method.

#### Parameters

- **node\_id** – the id of the node to write
- **value** – the value to write to the variable

```
class BaseCubeConfiguration(namespace_index: int = 3, polling_delay_sec: Union[int, float] = 5.0,  
polling_interval_sec: Union[int, float] = 1.0, timeout_status_change:  
Union[int, float] = 6, timeout_interval: Union[int, float] = 0.1,  
noise_level_measurement_channel_1: Union[int, float] = 100,  
noise_level_measurement_channel_2: Union[int, float] = 100,  
noise_level_measurement_channel_3: Union[int, float] = 100,  
noise_level_measurement_channel_4: Union[int, float] = 100)
```

Bases: object

Configuration dataclass for the BaseCube devices.

**clean\_values**()

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**namespace\_index: int = 3**

Namespace of the OPC variables, typically this is 3 (coming from Siemens)

**noise\_level\_measurement\_channel\_1: Union[int, float] = 100**

**noise\_level\_measurement\_channel\_2: Union[int, float] = 100**

**noise\_level\_measurement\_channel\_3: Union[int, float] = 100**

**noise\_level\_measurement\_channel\_4: Union[int, float] = 100**

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_delay\_sec:** `Union[int, float] = 5.0`

**polling\_interval\_sec:** `Union[int, float] = 1.0`

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**timeout\_interval:** `Union[int, float] = 0.1`

**timeout\_status\_change:** `Union[int, float] = 6`

**class BaseCubeOpcUaCommunication**(*config*)

Bases: `hvl_ccb.comm.opc.OpcUaCommunication`

Communication protocol specification for BaseCube devices.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class BaseCubeOpcUaCommunicationConfig**(*host: typing.Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], endpoint\_name: str = \_CubeOpcEndpoint.BASE\_CUBE, port: int = 4840, sub\_handler: hvl\_ccb.comm.opc.OpcUaSubHandler = <hvl\_ccb.dev.cube.base.\_BaseCubeSubscriptionHandler object>, update\_parameter: asyncua.ua.uaproTOCOL\_AUTO.CreateSubscriptionParameters = CreateSubscriptionParameters(RequestedPublishingInterval=1000, RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22, MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0), wait\_timeout\_retry\_sec: typing.Union[int, float] = 1, max\_timeout\_retry\_nr: int = 5)*)

Bases: `hvl_ccb.comm.opc.OpcUaCommunicationConfig`

Communication protocol configuration for OPC UA, specifications for the BaseCube devices.

**endpoint\_name:** `str = 'BaseCube'`

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**sub\_handler:** `hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.cube.base._BaseCubeSubscriptionHandler object>`

Subscription handler for data change events

### hvl\_ccb.dev.cube.constants

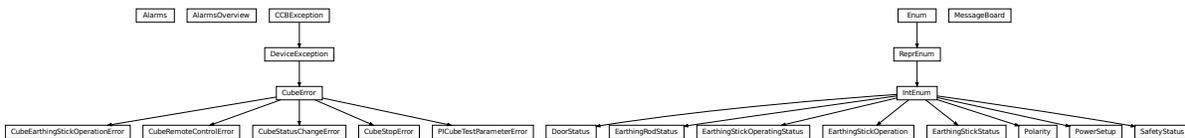

Constants, variable names for the BaseCube OPC-connected devices.

**class Alarms**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: `hvl_ccb.dev.cube.constants._AlarmEnumBase`

Alarms enumeration containing all variable NodeID strings for the alarm array.

**ALARM\_1** = "DB\_Alarm\_HMI"."Alarm1"

**ALARM\_10** = "DB\_Alarm\_HMI"."Alarm10"

**ALARM\_100** = "DB\_Alarm\_HMI"."Alarm100"

**ALARM\_101** = "DB\_Alarm\_HMI"."Alarm101"

**ALARM\_102** = "DB\_Alarm\_HMI"."Alarm102"

**ALARM\_103** = "DB\_Alarm\_HMI"."Alarm103"

**ALARM\_104** = "DB\_Alarm\_HMI"."Alarm104"

**ALARM\_105** = "DB\_Alarm\_HMI"."Alarm105"

**ALARM\_106** = "DB\_Alarm\_HMI"."Alarm106"

**ALARM\_107** = "DB\_Alarm\_HMI"."Alarm107"

**ALARM\_108** = "DB\_Alarm\_HMI"."Alarm108"

**ALARM\_109** = "DB\_Alarm\_HMI"."Alarm109"

**ALARM\_11** = "DB\_Alarm\_HMI"."Alarm11"

**ALARM\_110** = "DB\_Alarm\_HMI"."Alarm110"

**ALARM\_111** = "DB\_Alarm\_HMI"."Alarm111"

**ALARM\_112** = "DB\_Alarm\_HMI"."Alarm112"

**ALARM\_113** = "DB\_Alarm\_HMI"."Alarm113"

```
ALARM_114 = "DB_Alarm_HMI"."Alarm114"  
ALARM_115 = "DB_Alarm_HMI"."Alarm115"  
ALARM_116 = "DB_Alarm_HMI"."Alarm116"  
ALARM_117 = "DB_Alarm_HMI"."Alarm117"  
ALARM_118 = "DB_Alarm_HMI"."Alarm118"  
ALARM_119 = "DB_Alarm_HMI"."Alarm119"  
ALARM_12 = "DB_Alarm_HMI"."Alarm12"  
ALARM_120 = "DB_Alarm_HMI"."Alarm120"  
ALARM_121 = "DB_Alarm_HMI"."Alarm121"  
ALARM_122 = "DB_Alarm_HMI"."Alarm122"  
ALARM_123 = "DB_Alarm_HMI"."Alarm123"  
ALARM_124 = "DB_Alarm_HMI"."Alarm124"  
ALARM_125 = "DB_Alarm_HMI"."Alarm125"  
ALARM_126 = "DB_Alarm_HMI"."Alarm126"  
ALARM_127 = "DB_Alarm_HMI"."Alarm127"  
ALARM_128 = "DB_Alarm_HMI"."Alarm128"  
ALARM_129 = "DB_Alarm_HMI"."Alarm129"  
ALARM_13 = "DB_Alarm_HMI"."Alarm13"  
ALARM_130 = "DB_Alarm_HMI"."Alarm130"  
ALARM_131 = "DB_Alarm_HMI"."Alarm131"  
ALARM_132 = "DB_Alarm_HMI"."Alarm132"  
ALARM_133 = "DB_Alarm_HMI"."Alarm133"  
ALARM_134 = "DB_Alarm_HMI"."Alarm134"  
ALARM_135 = "DB_Alarm_HMI"."Alarm135"  
ALARM_136 = "DB_Alarm_HMI"."Alarm136"  
ALARM_137 = "DB_Alarm_HMI"."Alarm137"  
ALARM_138 = "DB_Alarm_HMI"."Alarm138"  
ALARM_139 = "DB_Alarm_HMI"."Alarm139"  
ALARM_14 = "DB_Alarm_HMI"."Alarm14"  
ALARM_140 = "DB_Alarm_HMI"."Alarm140"  
ALARM_141 = "DB_Alarm_HMI"."Alarm141"  
ALARM_142 = "DB_Alarm_HMI"."Alarm142"  
ALARM_143 = "DB_Alarm_HMI"."Alarm143"  
ALARM_144 = "DB_Alarm_HMI"."Alarm144"  
ALARM_145 = "DB_Alarm_HMI"."Alarm145"  
ALARM_146 = "DB_Alarm_HMI"."Alarm146"
```

```
ALARM_147 = 'DB_Alarm_HMI"."Alarm147"'
ALARM_148 = 'DB_Alarm_HMI"."Alarm148"'
ALARM_149 = 'DB_Alarm_HMI"."Alarm149"'
ALARM_15 = 'DB_Alarm_HMI"."Alarm15"'
ALARM_150 = 'DB_Alarm_HMI"."Alarm150"'
ALARM_151 = 'DB_Alarm_HMI"."Alarm151"'
ALARM_16 = 'DB_Alarm_HMI"."Alarm16"'
ALARM_17 = 'DB_Alarm_HMI"."Alarm17"'
ALARM_18 = 'DB_Alarm_HMI"."Alarm18"'
ALARM_19 = 'DB_Alarm_HMI"."Alarm19"'
ALARM_2 = 'DB_Alarm_HMI"."Alarm2"'
ALARM_20 = 'DB_Alarm_HMI"."Alarm20"'
ALARM_21 = 'DB_Alarm_HMI"."Alarm21"'
ALARM_22 = 'DB_Alarm_HMI"."Alarm22"'
ALARM_23 = 'DB_Alarm_HMI"."Alarm23"'
ALARM_24 = 'DB_Alarm_HMI"."Alarm24"'
ALARM_25 = 'DB_Alarm_HMI"."Alarm25"'
ALARM_26 = 'DB_Alarm_HMI"."Alarm26"'
ALARM_27 = 'DB_Alarm_HMI"."Alarm27"'
ALARM_28 = 'DB_Alarm_HMI"."Alarm28"'
ALARM_29 = 'DB_Alarm_HMI"."Alarm29"'
ALARM_3 = 'DB_Alarm_HMI"."Alarm3"'
ALARM_30 = 'DB_Alarm_HMI"."Alarm30"'
ALARM_31 = 'DB_Alarm_HMI"."Alarm31"'
ALARM_32 = 'DB_Alarm_HMI"."Alarm32"'
ALARM_33 = 'DB_Alarm_HMI"."Alarm33"'
ALARM_34 = 'DB_Alarm_HMI"."Alarm34"'
ALARM_35 = 'DB_Alarm_HMI"."Alarm35"'
ALARM_36 = 'DB_Alarm_HMI"."Alarm36"'
ALARM_37 = 'DB_Alarm_HMI"."Alarm37"'
ALARM_38 = 'DB_Alarm_HMI"."Alarm38"'
ALARM_39 = 'DB_Alarm_HMI"."Alarm39"'
ALARM_4 = 'DB_Alarm_HMI"."Alarm4"'
ALARM_40 = 'DB_Alarm_HMI"."Alarm40"'
ALARM_41 = 'DB_Alarm_HMI"."Alarm41"'
ALARM_42 = 'DB_Alarm_HMI"."Alarm42"'
```

```
ALARM_43 = '"DB_Alarm_HMI"."Alarm43"'
ALARM_44 = '"DB_Alarm_HMI"."Alarm44"'
ALARM_45 = '"DB_Alarm_HMI"."Alarm45"'
ALARM_46 = '"DB_Alarm_HMI"."Alarm46"'
ALARM_47 = '"DB_Alarm_HMI"."Alarm47"'
ALARM_48 = '"DB_Alarm_HMI"."Alarm48"'
ALARM_49 = '"DB_Alarm_HMI"."Alarm49"'
ALARM_5 = '"DB_Alarm_HMI"."Alarm5"'
ALARM_50 = '"DB_Alarm_HMI"."Alarm50"'
ALARM_51 = '"DB_Alarm_HMI"."Alarm51"'
ALARM_52 = '"DB_Alarm_HMI"."Alarm52"'
ALARM_53 = '"DB_Alarm_HMI"."Alarm53"'
ALARM_54 = '"DB_Alarm_HMI"."Alarm54"'
ALARM_55 = '"DB_Alarm_HMI"."Alarm55"'
ALARM_56 = '"DB_Alarm_HMI"."Alarm56"'
ALARM_57 = '"DB_Alarm_HMI"."Alarm57"'
ALARM_58 = '"DB_Alarm_HMI"."Alarm58"'
ALARM_59 = '"DB_Alarm_HMI"."Alarm59"'
ALARM_6 = '"DB_Alarm_HMI"."Alarm6"'
ALARM_60 = '"DB_Alarm_HMI"."Alarm60"'
ALARM_61 = '"DB_Alarm_HMI"."Alarm61"'
ALARM_62 = '"DB_Alarm_HMI"."Alarm62"'
ALARM_63 = '"DB_Alarm_HMI"."Alarm63"'
ALARM_64 = '"DB_Alarm_HMI"."Alarm64"'
ALARM_65 = '"DB_Alarm_HMI"."Alarm65"'
ALARM_66 = '"DB_Alarm_HMI"."Alarm66"'
ALARM_67 = '"DB_Alarm_HMI"."Alarm67"'
ALARM_68 = '"DB_Alarm_HMI"."Alarm68"'
ALARM_69 = '"DB_Alarm_HMI"."Alarm69"'
ALARM_7 = '"DB_Alarm_HMI"."Alarm7"'
ALARM_70 = '"DB_Alarm_HMI"."Alarm70"'
ALARM_71 = '"DB_Alarm_HMI"."Alarm71"'
ALARM_72 = '"DB_Alarm_HMI"."Alarm72"'
ALARM_73 = '"DB_Alarm_HMI"."Alarm73"'
ALARM_74 = '"DB_Alarm_HMI"."Alarm74"'
ALARM_75 = '"DB_Alarm_HMI"."Alarm75"'
```

```
ALARM_76 = '"DB_Alarm_HMI"."Alarm76"'
ALARM_77 = '"DB_Alarm_HMI"."Alarm77"'
ALARM_78 = '"DB_Alarm_HMI"."Alarm78"'
ALARM_79 = '"DB_Alarm_HMI"."Alarm79"'
ALARM_8 = '"DB_Alarm_HMI"."Alarm8"'
ALARM_80 = '"DB_Alarm_HMI"."Alarm80"'
ALARM_81 = '"DB_Alarm_HMI"."Alarm81"'
ALARM_82 = '"DB_Alarm_HMI"."Alarm82"'
ALARM_83 = '"DB_Alarm_HMI"."Alarm83"'
ALARM_84 = '"DB_Alarm_HMI"."Alarm84"'
ALARM_85 = '"DB_Alarm_HMI"."Alarm85"'
ALARM_86 = '"DB_Alarm_HMI"."Alarm86"'
ALARM_87 = '"DB_Alarm_HMI"."Alarm87"'
ALARM_88 = '"DB_Alarm_HMI"."Alarm88"'
ALARM_89 = '"DB_Alarm_HMI"."Alarm89"'
ALARM_9 = '"DB_Alarm_HMI"."Alarm9"'
ALARM_90 = '"DB_Alarm_HMI"."Alarm90"'
ALARM_91 = '"DB_Alarm_HMI"."Alarm91"'
ALARM_92 = '"DB_Alarm_HMI"."Alarm92"'
ALARM_93 = '"DB_Alarm_HMI"."Alarm93"'
ALARM_94 = '"DB_Alarm_HMI"."Alarm94"'
ALARM_95 = '"DB_Alarm_HMI"."Alarm95"'
ALARM_96 = '"DB_Alarm_HMI"."Alarm96"'
ALARM_97 = '"DB_Alarm_HMI"."Alarm97"'
ALARM_98 = '"DB_Alarm_HMI"."Alarm98"'
ALARM_99 = '"DB_Alarm_HMI"."Alarm99"'
```

```
class AlarmsOverview
```

```
    Bases: object
```

```
    Stores the status of all alarms / messages
```

```
    alarm_1 = 0
```

```
    alarm_10 = 0
```

```
    alarm_100 = 0
```

```
    alarm_101 = 0
```

```
    alarm_102 = 0
```

```
    alarm_103 = 0
```

```
    alarm_104 = 0
```

alarm\_105 = 0  
alarm\_106 = 0  
alarm\_107 = 0  
alarm\_108 = 0  
alarm\_109 = 0  
alarm\_11 = 0  
alarm\_110 = 0  
alarm\_111 = 0  
alarm\_112 = 0  
alarm\_113 = 0  
alarm\_114 = 0  
alarm\_115 = 0  
alarm\_116 = 0  
alarm\_117 = 0  
alarm\_118 = 0  
alarm\_119 = 0  
alarm\_12 = 0  
alarm\_120 = 0  
alarm\_121 = 0  
alarm\_122 = 0  
alarm\_123 = 0  
alarm\_124 = 0  
alarm\_125 = 0  
alarm\_126 = 0  
alarm\_127 = 0  
alarm\_128 = 0  
alarm\_129 = 0  
alarm\_13 = 0  
alarm\_130 = 0  
alarm\_131 = 0  
alarm\_132 = 0  
alarm\_133 = 0  
alarm\_134 = 0  
alarm\_135 = 0  
alarm\_136 = 0  
alarm\_137 = 0

alarm\_138 = 0  
alarm\_139 = 0  
alarm\_14 = 0  
alarm\_140 = 0  
alarm\_141 = 0  
alarm\_142 = 0  
alarm\_143 = 0  
alarm\_144 = 0  
alarm\_145 = 0  
alarm\_146 = 0  
alarm\_147 = 0  
alarm\_148 = 0  
alarm\_149 = 0  
alarm\_15 = 0  
alarm\_150 = 0  
alarm\_151 = 0  
alarm\_16 = 0  
alarm\_17 = 0  
alarm\_18 = 0  
alarm\_19 = 0  
alarm\_2 = 0  
alarm\_20 = 0  
alarm\_21 = 0  
alarm\_22 = 0  
alarm\_23 = 0  
alarm\_24 = 0  
alarm\_25 = 0  
alarm\_26 = 0  
alarm\_27 = 0  
alarm\_28 = 0  
alarm\_29 = 0  
alarm\_3 = 0  
alarm\_30 = 0  
alarm\_31 = 0  
alarm\_32 = 0  
alarm\_33 = 0

alarm\_34 = 0  
alarm\_35 = 0  
alarm\_36 = 0  
alarm\_37 = 0  
alarm\_38 = 0  
alarm\_39 = 0  
alarm\_4 = 0  
alarm\_40 = 0  
alarm\_41 = 0  
alarm\_42 = 0  
alarm\_43 = 0  
alarm\_44 = 0  
alarm\_45 = 0  
alarm\_46 = 0  
alarm\_47 = 0  
alarm\_48 = 0  
alarm\_49 = 0  
alarm\_5 = 0  
alarm\_50 = 0  
alarm\_51 = 0  
alarm\_52 = 0  
alarm\_53 = 0  
alarm\_54 = 0  
alarm\_55 = 0  
alarm\_56 = 0  
alarm\_57 = 0  
alarm\_58 = 0  
alarm\_59 = 0  
alarm\_6 = 0  
alarm\_60 = 0  
alarm\_61 = 0  
alarm\_62 = 0  
alarm\_63 = 0  
alarm\_64 = 0  
alarm\_65 = 0  
alarm\_66 = 0

alarm\_67 = 0  
alarm\_68 = 0  
alarm\_69 = 0  
alarm\_7 = 0  
alarm\_70 = 0  
alarm\_71 = 0  
alarm\_72 = 0  
alarm\_73 = 0  
alarm\_74 = 0  
alarm\_75 = 0  
alarm\_76 = 0  
alarm\_77 = 0  
alarm\_78 = 0  
alarm\_79 = 0  
alarm\_8 = 0  
alarm\_80 = 0  
alarm\_81 = 0  
alarm\_82 = 0  
alarm\_83 = 0  
alarm\_84 = 0  
alarm\_85 = 0  
alarm\_86 = 0  
alarm\_87 = 0  
alarm\_88 = 0  
alarm\_89 = 0  
alarm\_9 = 0  
alarm\_90 = 0  
alarm\_91 = 0  
alarm\_92 = 0  
alarm\_93 = 0  
alarm\_94 = 0  
alarm\_95 = 0  
alarm\_96 = 0  
alarm\_97 = 0  
alarm\_98 = 0  
alarm\_99 = 0

**exception CubeEarthingStickOperationError**

Bases: *hvl\_ccb.dev.cube.constants.CubeError*

**exception CubeError**

Bases: *hvl\_ccb.dev.base.DeviceException*

**exception CubeRemoteControlError**

Bases: *hvl\_ccb.dev.cube.constants.CubeError*

**exception CubeStatusChangeError**

Bases: *hvl\_ccb.dev.cube.constants.CubeError*

**exception CubeStopError**

Bases: *hvl\_ccb.dev.cube.constants.CubeError*

**class DoorStatus**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *aenum.IntEnum*

Possible status values for doors.

**CLOSED = 2**

Door is closed, but not locked.

**ERROR = 4**

Door has an error or was opened in locked state (either with emergency stop or from the inside).

**INACTIVE = 0**

not enabled in BaseCube HMI setup, this door is not supervised.

**LOCKED = 3**

Door is closed and locked (safe state).

**OPEN = 1**

Door is open.

**class EarthingRodStatus**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *aenum.IntEnum*

Possible status values for earthing rods.

**EXPERIMENT\_BLOCKED = 0**

earthing rod is somewhere in the experiment and blocks the start of the experiment

**EXPERIMENT\_READY = 1**

earthing rod is hanging next to the door, experiment is ready to operate

**class EarthingStickOperatingStatus**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *aenum.IntEnum*

Operating Status for an earthing stick. Stick can be used in auto or manual mode.

**AUTO = 0**

**MANUAL = 1**

**class EarthingStickOperation**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *aenum.IntEnum*

Operation of the earthing stick in manual operating mode. Can be closed or opened.

**CLOSE = 1**

**OPEN = 0**

**class EarthingStickStatus**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*,  
*boundary=None*)

Bases: `aenum.IntEnum`

Status of an earthing stick. These are the possible values in the status integer e.g. in `_EarthingStick.status`.

**CLOSED = 1**

**ERROR = 3**

**INACTIVE = 0**

**OPEN = 2**

**class MessageBoard**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.dev.cube.constants._LineEnumBase`

Variable NodeID strings for message board lines.

**LINE\_1 = "DB\_OPC\_Connection"."Is\_status\_Line\_1"**

**LINE\_10 = "DB\_OPC\_Connection"."Is\_status\_Line\_10"**

**LINE\_11 = "DB\_OPC\_Connection"."Is\_status\_Line\_11"**

**LINE\_12 = "DB\_OPC\_Connection"."Is\_status\_Line\_12"**

**LINE\_13 = "DB\_OPC\_Connection"."Is\_status\_Line\_13"**

**LINE\_14 = "DB\_OPC\_Connection"."Is\_status\_Line\_14"**

**LINE\_15 = "DB\_OPC\_Connection"."Is\_status\_Line\_15"**

**LINE\_2 = "DB\_OPC\_Connection"."Is\_status\_Line\_2"**

**LINE\_3 = "DB\_OPC\_Connection"."Is\_status\_Line\_3"**

**LINE\_4 = "DB\_OPC\_Connection"."Is\_status\_Line\_4"**

**LINE\_5 = "DB\_OPC\_Connection"."Is\_status\_Line\_5"**

**LINE\_6 = "DB\_OPC\_Connection"."Is\_status\_Line\_6"**

**LINE\_7 = "DB\_OPC\_Connection"."Is\_status\_Line\_7"**

**LINE\_8 = "DB\_OPC\_Connection"."Is\_status\_Line\_8"**

**LINE\_9 = "DB\_OPC\_Connection"."Is\_status\_Line\_9"**

**exception PICubeTestParameterError**

Bases: `hvl_ccb.dev.cube.constants.CubeError`

**class Polarity**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `aenum.IntEnum`

An enumeration.

**NEGATIVE = 0**

**POSITIVE = 1**

**class PowerSetup**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `aenum.IntEnum`

Possible power setups corresponding to the value of variable `Power.setup`. The values for `slope_min` are experimentally defined, below these values the slope is more like a staircase

The name of the first argument needs to be 'value', otherwise the IntEnum is not working correctly.

AC\_100KV = 3

AC\_150KV = 4

AC\_200KV = 5

AC\_50KV = 2

DC\_140KV = 7

DC\_280KV = 8

EXTERNAL\_SOURCE = 1

IMPULSE\_140KV = 9

NO\_SOURCE = 0

POWER\_INVERTER\_220V = 6

STOP\_SAFETY\_STATUSES: Tuple[[hvl\\_ccb.dev.cube.constants.SafetyStatus](#), ...] =  
(<SafetyStatus.GREEN\_NOT\_READY: 1>, <SafetyStatus.GREEN\_READY: 2>)

BaseCube's safety statuses required to close the connection to the device.

**class SafetyStatus**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: aenum.IntEnum

Safety status values that are possible states returned from [hvl\\_ccb.dev.cube.base.BaseCube.status\(\)](#).  
These values correspond to the states of the BaseCube's safety circuit statemachine.

ERROR = 6

GREEN\_NOT\_READY = 1

GREEN\_READY = 2


INITIALIZING = 0

QUICK\_STOP = 5

RED\_OPERATE = 4

RED\_READY = 3

## hvl\_ccb.dev.cube.picube

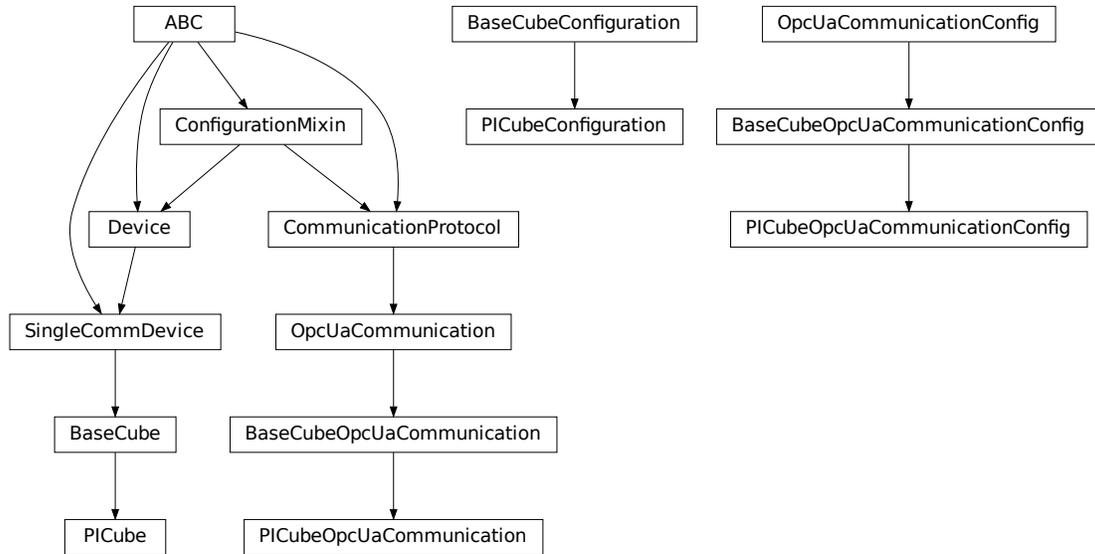

A PICube is a BaseCube with build in Power Inverter

**class** `PICube`(*com*, *dev\_config=None*)

Bases: `hvl_ccb.dev.cube.base.BaseCube`

Variant of the BaseCube with build in Power Inverter

**static** `config_cls()`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property** `current_primary: float`

Read the current primary current at the output of the frequency converter (before transformer).

**Returns** primary current in A

**static** `default_com_cls()`

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**property** `frequency: float`

Read the electrical frequency of the current PICube setup.

**Returns** the frequency in Hz

**property** `operate: Optional[bool]`

Indicates if 'operate' is activated. 'operate' means locket safety circuit, red lamps, high voltage on and locked safety switches.

**Returns** *True* if operate is activated (RedOperate), *False* if ready is deactivated (RedReady), *None* otherwise

**property** `polarity: Optional[hvl_ccb.dev.cube.constants.Polarity]`

Polarity of a DC setup. :return: if a DC setup is programmed the polarity is returned, else None.

**property power\_setup:** [hvl\\_ccb.dev.cube.constants.PowerSetup](#)

Return the power setup selected in the PICube's settings.

**Returns** the power setup

**property voltage\_actual:** **float**

Reads the actual measured voltage and returns the value in V.

**Returns** the actual voltage of the setup in V.

**property voltage\_max:** **float**

Reads the maximum voltage of the setup and returns in V.

**Returns** the maximum voltage of the setup in V.

**property voltage\_primary:** **float**

Read the current primary voltage at the output of the frequency converter (before transformer).

**Returns** primary voltage in V

```
class PICubeConfiguration(namespace_index: int = 3, polling_delay_sec: Union[int, float] = 5.0,
                          polling_interval_sec: Union[int, float] = 1.0, timeout_status_change: Union[int,
                          float] = 6, timeout_interval: Union[int, float] = 0.1,
                          noise_level_measurement_channel_1: Union[int, float] = 100,
                          noise_level_measurement_channel_2: Union[int, float] = 100,
                          noise_level_measurement_channel_3: Union[int, float] = 100,
                          noise_level_measurement_channel_4: Union[int, float] = 100,
                          timeout_test_parameters: 'Number' = 2.0)
```

Bases: [hvl\\_ccb.dev.cube.base.BaseCubeConfiguration](#)

**clean\_values()**

**force\_value**(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**timeout\_test\_parameters:** **Union[int, float] = 2.0**

```
class PICubeOpcUaCommunication(config)
```

Bases: [hvl\\_ccb.dev.cube.base.BaseCubeOpcUaCommunication](#)

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class PICubeOpcUaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],  
endpoint_name: 'str' = <_CubeOpcEndpoint.PI_CUBE: 'PICube'>,  
port: int = 4840, sub_handler:  
hvl_ccb.comm.opc.OpcUaSubHandler =  
<hvl_ccb.dev.cube.base._BaseCubeSubscriptionHandler object at  
0x7f7a2b565750>, update_parameter:  
asynqua.ua.uaprotoocol_auto.CreateSubscriptionParameters =  
CreateSubscriptionParameters(RequestedPublishingInterval=1000,  
RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22,  
MaxNotificationsPerPublish=10000, PublishingEnabled=True,  
Priority=0), wait_timeout_retry_sec: Union[int, float] = 1,  
max_timeout_retry_nr: int = 5)
```

Bases: [hvl\\_ccb.dev.cube.base.BaseCubeOpcUaCommunicationConfig](#)

**endpoint\_name: str = 'PICube'**

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


## Module contents

Cube package with implementation for system versions from 2019 on (new concept with hard-PLC Siemens S7-1500 as CPU).

### hvl\_ccb.dev.highland\_t560

#### Submodules

#### hvl\_ccb.dev.highland\_t560.base

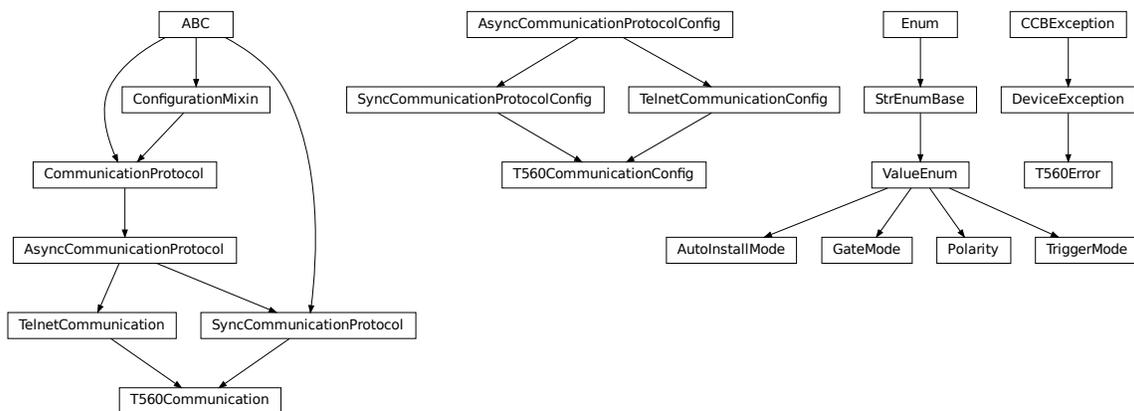

Module containing base device and communication classes and enums.

Communication with device is performed via its ethernet port and a Telnet connection.

**class AutoInstallMode**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Modes for installing configuration settings to the device.

**INSTALL = 1**

**OFF = 0**

**QUEUE = 2**

**class GateMode**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Available T560 gate modes

**INPUT = 'INP'**

**OFF = 'OFF'**

**OUTPUT = 'OUT'**

**class Polarity**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Possible channel polarity states

**ACTIVE\_HIGH** = 'POS'

**ACTIVE\_LOW** = 'NEG'

**class T560Communication**(*configuration*)

Bases: [hvl\\_ccb.comm.base.SyncCommunicationProtocol](#), [hvl\\_ccb.comm.telnet.TelnetCommunication](#)

Communication class for T560. It uses a TelnetCommunication with the SyncCommunicationProtocol

**static config\_cls**()

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query**(*command: str*) → str

Send a command to the device and handle the response.

For device setting queries, response will be 'OK' if successful, or '??' if setting cannot be carried out, raising an error.

**Parameters** **command** – Command string to be sent

**Raises** [T560Error](#) – if no response is received, or if the device responds with an error message.

**Returns** Response from the device.

**class T560CommunicationConfig**(*terminator: bytes = b'\r', encoding: str = 'utf-8', encoding\_error\_handling: str = 'strict', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10, host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address, NoneType] = None, port: int = 2000, timeout: Union[int, float] = 0.2)*)

Bases: [hvl\\_ccb.comm.base.SyncCommunicationProtocolConfig](#), [hvl\\_ccb.comm.telnet.TelnetCommunicationConfig](#)

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 2000**

Port at which the host is listening

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**terminator:** `bytes = b'\r'`

The terminator character. Typically this is `b'\r\n'` or `b'\n'`, but can also be `b'\r'` or other combinations. This defines the end of a single line.

**exception** `T560Error`

Bases: `hvl_ccb.dev.base.DeviceException`

T560 related errors.

**class** `TriggerMode`(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.ValueEnum`

Available T560 trigger modes

`COMMAND = 'REM'`

`EXT_FALLING_EDGE = 'NEG'`

`EXT_RISING_EDGE = 'POS'`


`INT_SYNTHESIZER = 'SYN'`

`OFF = 'OFF'`

**`hvl_ccb.dev.highland_t560.channel`**

Module for controlling pulse output channels A, B, C and D.

## hvl\_ccb.dev.highland\_t560.device

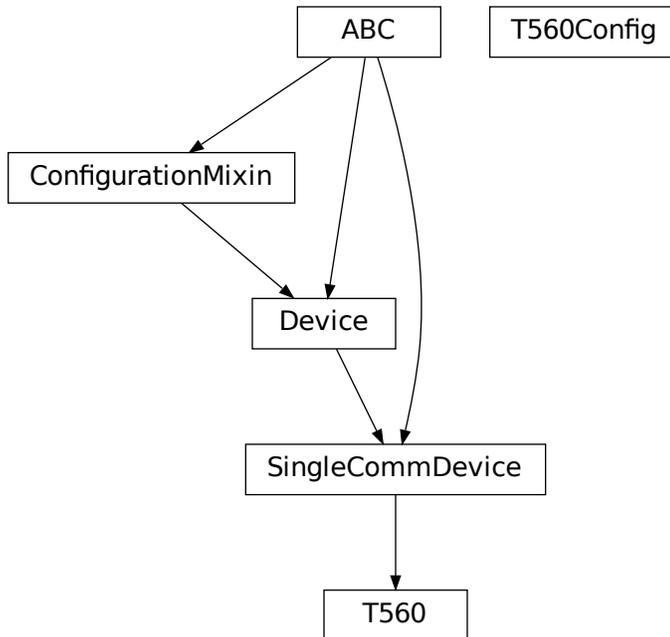

including TRIG, CLOCK and GATE I/Os.

Module for controlling device,

**class T560**(*com, dev\_config=None*)

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

**activate\_clock\_output**()

Outputs 10 MHz clock signal

**property auto\_install\_mode:** *hvl\_ccb.dev.highland\_t560.base.AutoInstallMode*

Check the autoinstall settings of the T560. The autoinstall mode sets how changes to device settings are applied. See manual section 4.7.2 for more information about these modes.

**property ch\_a:** *hvl\_ccb.dev.highland\_t560.channel.\_Channel*

Channel A of T560

**property ch\_b:** *hvl\_ccb.dev.highland\_t560.channel.\_Channel*

Channel B of T560

**property ch\_c:** *hvl\_ccb.dev.highland\_t560.channel.\_Channel*

Channel C of T560

**property ch\_d:** *hvl\_ccb.dev.highland\_t560.channel.\_Channel*

Channel D of T560

**static config\_cls**()

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**disarm\_trigger()**

Disarm DDG by disabling all trigger sources.

**fire\_trigger()**

Fire a software trigger.

**property frequency: float**

The frequency of the timing cycle in Hz.

**property gate\_mode: *hvl\_ccb.dev.highland\_t560.base.GateMode***

Check the mode setting of the GATE I/O port.

**property gate\_polarity: *hvl\_ccb.dev.highland\_t560.base.Polarity***

Check the polarity setting of the GATE I/O port.

**load\_device\_configuration()**

Load the settings saved in nonvolatile memory.

**property period: float**

The period of the timing cycle (time between triggers) in seconds.

**save\_device\_configuration()**

Save the current settings to nonvolatile memory.

**property trigger\_level**

Get external trigger level.

**property trigger\_mode**

Get device trigger source.

**use\_external\_clock()**

Finds and accepts an external clock signal to the CLOCK input

**class T560Config**

Bases: object

**auto\_install\_mode = 1**

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

## Module contents

This module establishes methods for interfacing with the Highland Technology T560-2 via its ethernet adapter with a telnet communication protocol.

The T560 is a small digital delay & pulse generator. It outputs up to four individually timed pulses with 10-ps precision, given an internal or external trigger.


This module introduces methods for configuring channels, gating, and triggering. Further documentation and a more extensive command list may be obtained from:

<https://www.highlandtechnology.com/DSS/T560DS.shtml>

## hvl\_ccb.dev.supercube2015

### Submodules

#### hvl\_ccb.dev.supercube2015.base

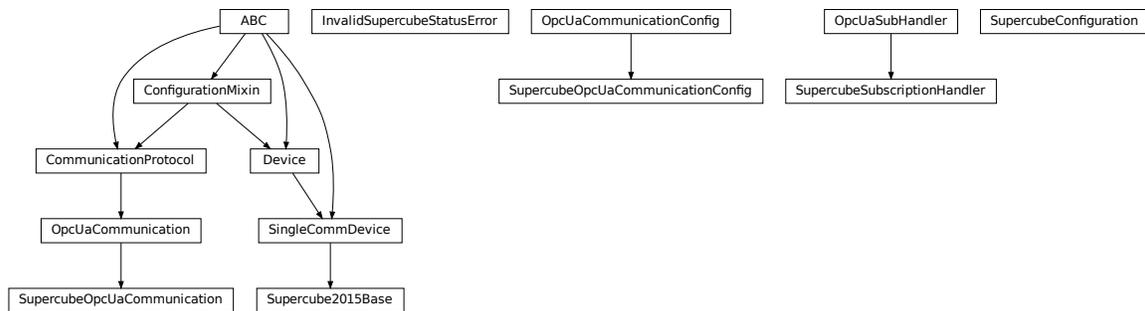

Base classes for the Supercube device.

**exception InvalidSupercubeStatusError**

Bases: Exception

Exception raised when supercube has invalid status.

**class Supercube2015Base**(com, dev\_config=None)

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Base class for Supercube variants.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**get\_cee16\_socket()** → bool

Read the on-state of the IEC CEE16 three-phase power socket.

**Returns** the on-state of the CEE16 power socket

**get\_door\_status(door: int)** → *hvl\_ccb.dev.supercube2015.constants.DoorStatus*

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

**Parameters** `door` – the door number (1..3)

**Returns** the door status

**get\_earthing\_manual(number: int)** → bool

Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** earthing stick manual status

**get\_earthing\_status(number: int)** → int

Get the status of an earthing stick, whether it is closed, open or undefined (moving).

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** earthing stick status; see `constants.EarthingStickStatus`

**get\_measurement\_ratio(channel: int)** → float

Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

**Attention:** Supercube 2015 does not have a separate ratio for every analog input. Therefore there is only one ratio for `channel = 1`.

**Parameters** `channel` – number of the input channel (1..4)

**Returns** the ratio

**get\_measurement\_voltage(channel: int)** → float

Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

**Attention:** In contrast to the *new* Supercube, the old one returns here the input voltage read at the ADC. It is not scaled by a factor.

**Parameters** `channel` – number of the input channel (1..4)

**Returns** measured voltage

**get\_status()** → int

Get the safety circuit status of the Supercube.

**Returns** the safety status of the supercube's state machine; see `constants.SafetyStatus`.

**get\_support\_input(port: int, contact: int)** → bool

Get the state of a support socket input.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

**Returns** digital input read state

**get\_support\_output**(*port: int, contact: int*) → bool  
Get the state of a support socket output.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

**Returns** digital output read state

**get\_t13\_socket**(*port: int*) → bool  
Read the state of a SEV T13 power socket.

**Parameters** **port** – is the socket number, one of *constants.T13\_SOCKET\_PORTS*

**Returns** on-state of the power socket

**horn**(*state: bool*) → *None*  
Turns acoustic horn on or off.

**Parameters** **state** – Turns horn on (True) or off (False)

**operate**(*state: bool*) → *None*  
Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

**Parameters** **state** – set operate state

**quit\_error**() → *None*  
Quits errors that are active on the Supercube.

**read**(*node\_id: str*)  
Local wrapper for the OPC UA communication protocol read method.

**Parameters** **node\_id** – the id of the node to read.

**Returns** the value of the variable

**ready**(*state: bool*) → *None*  
Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

**Parameters** **state** – set ready state

**set\_cee16\_socket**(*state: bool*) → *None*  
Switch the IEC CEE16 three-phase power socket on or off.

**Parameters** **state** – desired on-state of the power socket

**Raises** **TypeError** – if state is not of type bool

**set\_earthing\_manual**(*number: int, manual: bool*) → *None*  
Set the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

**Parameters**

- **number** – number of the earthing stick (1..6)
- **manual** – earthing stick manual status (True or False)

**set\_remote\_control**(*state: bool*) → *None*

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

**Parameters** **state** – desired remote control state

**set\_support\_output**(*port: int, contact: int, state: bool*) → *None*

Set the state of a support output socket.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **state** – is the desired state of the support output

**Raises** **TypeError** – when state is not of type bool

**set\_support\_output\_impulse**(*port: int, contact: int, duration: float = 0.2, pos\_pulse: bool = True*) → *None*

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the *pos\_pulse* argument.

This function is blocking.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **duration** – is the length of the impulse in seconds
- **pos\_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

**Raises** **TypeError** – when state is not of type bool

**set\_t13\_socket**(*port: int, state: bool*) → *None*

Set the state of a SEV T13 power socket.

**Parameters**

- **port** – is the socket number, one of *constants.T13\_SOCKET\_PORTS*
- **state** – is the desired on-state of the socket

**Raises**

- **ValueError** – when port is not valid
- **TypeError** – when state is not of type bool

**start**() → *None*

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

**stop**() → *None*

Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

**write**(*node\_id, value*) → *None*

Local wrapper for the OPC UA communication protocol write method.

**Parameters**

- **node\_id** – the id of the node to read
- **value** – the value to write to the variable

**class SupercubeConfiguration**(*namespace\_index: int = 7*)

Bases: object

Configuration dataclass for the Supercube devices.

**clean\_values**()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**namespace\_index: int = 7**

Namespace of the OPC variables, typically this is 3 (coming from Siemens)

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**class SupercubeOpcUaCommunication**(*config*)

Bases: *hvl\_ccb.comm.opc.OpcUaCommunication*

Communication protocol specification for Supercube devices.

**static config\_cls**()

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class SupercubeOpcUaCommunicationConfig(host: typing.Union[str, ipaddress.IPv4Address,
                                         ipaddress.IPv6Address], endpoint_name: str, port: int = 4845,
                                         sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
                                         <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler
                                         object>, update_parameter:
                                         asyncua.ua.uaprotoool_auto.CreateSubscriptionParameters =
                                         CreateSubscriptionParameters(
                                         RequestedPublishingInterval=1000,
                                         RequestedLifetimeCount=300,
                                         RequestedMaxKeepAliveCount=22,
                                         MaxNotificationsPerPublish=10000, PublishingEnabled=True,
                                         Priority=0), wait_timeout_retry_sec: typing.Union[int, float] =
                                         1, max_timeout_retry_nr: int = 5)
```

Bases: `hvl_ccb.comm.opc.OpcUaCommunicationConfig`

Communication protocol configuration for OPC UA, specifications for the Supercube devices.

**force\_value**(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 4845**

Port of the OPC UA server to connect to.

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**sub\_handler: hvl\_ccb.comm.opc.OpcUaSubHandler =**  
**<hvl\_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object>**

Subscription handler for data change events

**class SupercubeSubscriptionHandler**

Bases: `hvl_ccb.comm.opc.OpcUaSubHandler`

OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.


**datachange\_notification**(node: asyncua.sync.SyncNode, val, data)

In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.

### Parameters

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

### hvl\_ccb.dev.supercube2015.constants

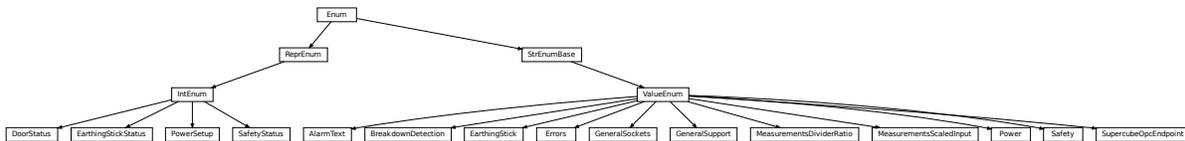

Constants, variable names for the Supercube OPC-connected devices.

**class AlarmText** (*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)  
 Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the [AlarmText.get\(\)](#) method to retrieve the enum of an alarm number.

```

Alarm0 = 'No Alarm.'
Alarm1 = 'STOP Safety switch 1 error'
Alarm10 = 'STOP Earthing stick 2 error'
Alarm11 = 'STOP Earthing stick 3 error'
Alarm12 = 'STOP Earthing stick 4 error'
Alarm13 = 'STOP Earthing stick 5 error'
Alarm14 = 'STOP Earthing stick 6 error'
Alarm17 = 'STOP Source switch error'
Alarm19 = 'STOP Fence 1 error'
Alarm2 = 'STOP Safety switch 2 error'
Alarm20 = 'STOP Fence 2 error'
Alarm21 = 'STOP Control error'
Alarm22 = 'STOP Power outage'
Alarm3 = 'STOP Emergency Stop 1'
Alarm4 = 'STOP Emergency Stop 2'
Alarm5 = 'STOP Emergency Stop 3'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error'
  
```

**classmethod** `get(alarm: int)`

Get the attribute of this enum for an alarm number.

**Parameters** `alarm` – the alarm number

**Returns** the enum for the desired alarm number

`not_defined = 'NO ALARM TEXT DEFINED'`

**class** `BreakdownDetection(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)`

Bases: `hvl_ccb.utils.enum.ValueEnum`

Node ID strings for the breakdown detection.

**activated = 'hvl-ipc.WINAC.SYSTEM\_COMPONENTS.Breakdowndetection.connect'**

Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the system or not.

**reset = 'hvl-ipc.WINAC.Support60uTA'**

Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

**triggered = 'hvl-ipc.WINAC.SYSTEM\_COMPONENTS.Breakdowndetection.triggered'**

Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using the safety circuit state, therefore no method is implemented to read this out directly.

**class** `DoorStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)`

Bases: `aenum.IntEnum`

Possible status values for doors.

**closed = 2**

Door is closed, but not locked.

**error = 4**

Door has an error or was opened in locked state (either with emergency stop or from the inside).

**inactive = 0**

not enabled in Supercube HMI setup, this door is not supervised.

**locked = 3**

Door is closed and locked (safe state).

**open = 1**

Door is open.

**class** `EarthingStick(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)`

Bases: `hvl_ccb.utils.enum.ValueEnum`

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the earthing in manual mode.

**classmethod** `manual(number: int)`

Get the manual enum attribute for an earthing stick number.

**Parameters** `number` – the earthing stick (1..6)

**Returns** the manual enum

`manual_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.MANUAL'`

`manual_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.MANUAL'`

`manual_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.MANUAL'`

`manual_4 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.MANUAL'`

```
manual_5 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.MANUAL'  
manual_6 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.MANUAL'  
status_1_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.CLOSE'  
status_1_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.CONNECT'  
status_1_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.OPEN'  
status_2_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.CLOSE'  
status_2_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.CONNECT'  
status_2_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.OPEN'  
status_3_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.CLOSE'  
status_3_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.CONNECT'  
status_3_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.OPEN'  
status_4_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.CLOSE'  
status_4_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.CONNECT'  
status_4_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.OPEN'  
status_5_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.CLOSE'  
status_5_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.CONNECT'  
status_5_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.OPEN'  
status_6_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.CLOSE'  
status_6_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.CONNECT'  
status_6_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.OPEN'
```

**classmethod** `status_closed(number: int)`

Get the status enum attribute for an earthing stick number.

**Parameters** `number` – the earthing stick (1..6)

**Returns** the status enum

**classmethod** `status_connected(number: int)`

Get the status enum attribute for an earthing stick number.

**Parameters** `number` – the earthing stick (1..6)

**Returns** the status enum

**classmethod** `status_open(number: int)`

Get the status enum attribute for an earthing stick number.

**Parameters** `number` – the earthing stick (1..6)

**Returns** the status enum

**class** `EarthingStickStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)`

Bases: `aenum.IntEnum`

Status of an earthing stick. These are the possible values in the status integer e.g. in `EarthingStick.status_1`.

**closed = 1**

Earthing is closed (safe).

**error = 3**

Earthing is in error, e.g. when the stick did not close correctly or could not open.

**inactive = 0**

Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be enabled in the Supercube HMI setup.

**open = 2**

Earthing is open (not safe).

**class Errors**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Variable NodeID strings for information regarding error, warning and message handling.

**quit = 'hvl-ipc.WINAC.SYSTEMSTATE.Faultconfirmation'**

Writable boolean for the error quit button.

**stop = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'**

Boolean read-only variable telling if a stop is active.

**stop\_number = 'hvl-ipc.WINAC.SYSTEMSTATE.Errornumber'**

**class GeneralSockets**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

NodeID strings for the power sockets (3x T13 and 1xCEE16).

**cee16 = 'hvl-ipc.WINAC.SYSTEM\_COMPONENTS.CEE16'**

CEE16 socket (writeable boolean).

**t13\_1 = 'hvl-ipc.WINAC.SYSTEM\_COMPONENTS.T13\_1'**

SEV T13 socket No. 1 (writeable boolean).

**t13\_2 = 'hvl-ipc.WINAC.SYSTEM\_COMPONENTS.T13\_2'**

SEV T13 socket No. 2 (writeable boolean).

**t13\_3 = 'hvl-ipc.WINAC.SYSTEM\_COMPONENTS.T13\_3'**

SEV T13 socket No. 3 (writeable boolean).

**class GeneralSupport**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

NodeID strings for the support inputs and outputs.

**in\_1\_1 = 'hvl-ipc.WINAC.Support1InA'**

**in\_1\_2 = 'hvl-ipc.WINAC.Support1InB'**

**in\_2\_1 = 'hvl-ipc.WINAC.Support2InA'**

**in\_2\_2 = 'hvl-ipc.WINAC.Support2InB'**

**in\_3\_1 = 'hvl-ipc.WINAC.Support3InA'**

**in\_3\_2 = 'hvl-ipc.WINAC.Support3InB'**

**in\_4\_1 = 'hvl-ipc.WINAC.Support4InA'**

**in\_4\_2 = 'hvl-ipc.WINAC.Support4InB'**

**in\_5\_1 = 'hvl-ipc.WINAC.Support5InA'**

**in\_5\_2 = 'hvl-ipc.WINAC.Support5InB'**

```
in_6_1 = 'hvl-ipc.WINAC.Support6InA'
```

```
in_6_2 = 'hvl-ipc.WINAC.Support6InB'
```

```
classmethod input(port, contact)
```

Get the NodeID string for a support input.

**Parameters**

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

**Returns** the node id string

```
out_1_1 = 'hvl-ipc.WINAC.Support10OutA'
```

```
out_1_2 = 'hvl-ipc.WINAC.Support10OutB'
```

```
out_2_1 = 'hvl-ipc.WINAC.Support20OutA'
```

```
out_2_2 = 'hvl-ipc.WINAC.Support20OutB'
```

```
out_3_1 = 'hvl-ipc.WINAC.Support30OutA'
```

```
out_3_2 = 'hvl-ipc.WINAC.Support30OutB'
```

```
out_4_1 = 'hvl-ipc.WINAC.Support40OutA'
```

```
out_4_2 = 'hvl-ipc.WINAC.Support40OutB'
```

```
out_5_1 = 'hvl-ipc.WINAC.Support50OutA'
```

```
out_5_2 = 'hvl-ipc.WINAC.Support50OutB'
```

```
out_6_1 = 'hvl-ipc.WINAC.Support60OutA'
```

```
out_6_2 = 'hvl-ipc.WINAC.Support60OutB'
```

```
classmethod output(port, contact)
```

Get the NodeID string for a support output.

**Parameters**

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

**Returns** the node id string

```
class MeasurementsDividerRatio(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

```
classmethod get(channel: int)
```

Get the attribute for an input number.

**Parameters** **channel** – the channel number (1..4)

**Returns** the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.DivididerRatio'
```

```
class MeasurementsScaledInput(value=<no_arg>, names=None, module=None, type=None, start=1,
                               boundary=None)
```

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in [MeasurementsDividerRatio](#).

```
classmethod get(channel: int)
```

Get the attribute for an input number.

**Parameters** **channel** – the channel number (1..4)

**Returns** the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI1Volt'
```

```
input_2 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI2Volt'
```

```
input_3 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI3Volt'
```

```
input_4 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI4Volt'
```

```
class Power(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Variable NodeID strings concerning power data.

```
current_primary = 'hvl-ipc.WINAC.SYSTEM_INTERN.FUCurrentprim'
```

Primary current in ampere, measured by the frequency converter. (read-only)

```
frequency = 'hvl-ipc.WINAC.FU.Frequency'
```

Frequency converter output frequency. (read-only)

```
setup = 'hvl-ipc.WINAC.FU.TrafoSetup'
```

Power setup that is configured using the Supercube HMI. The value corresponds to the ones in [PowerSetup](#). (read-only)

```
voltage_max = 'hvl-ipc.WINAC.FU.maxVoltagekV'
```

Maximum voltage allowed by the current experimental setup. (read-only)

```
voltage_primary = 'hvl-ipc.WINAC.SYSTEM_INTERN.FUVoltageprim'
```

Primary voltage in volts, measured by the frequency converter at its output. (read-only)

```
voltage_slope = 'hvl-ipc.WINAC.FU.dUdt_-1'
```

Voltage slope in V/s.

```
voltage_target = 'hvl-ipc.WINAC.FU.SOLL'
```

Target voltage setpoint in V.

```
class PowerSetup(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: [aenum.IntEnum](#)

Possible power setups corresponding to the value of variable [Power.setup](#).

```
AC_DoubleStage_150kV = 3
```

AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum voltage of 150kV.

```
AC_DoubleStage_200kV = 4
```

AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

```
AC_SingleStage_100kV = 2
```

AC voltage with MWB transformer set to 100kV maximum voltage.

**AC\_SingleStage\_50kV = 1**

AC voltage with MWB transformer set to 50kV maximum voltage.

**DC\_DoubleStage\_280kV = 7**

DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single stage transformer with Greinacher voltage doubling rectifier)

**DC\_SingleStage\_140kV = 6**

DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

**External = 0**

External power supply fed through blue CEE32 input using isolation transformer and safety switches of the Supercube, or using an external safety switch attached to the Supercube Type B.

**Internal = 5**

Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

**class Safety**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

**horn = 'hvl-ipc.WINAC.SYSTEM\_INTERN.hornen'**

Writable boolean to manually turn on or off the horn

**status\_error = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'**

**status\_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN'**

**status\_ready\_for\_red = 'hvl-ipc.WINAC.SYSTEMSTATE.ReadyForRed'**

Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in [SafetyStatus](#).

**status\_red = 'hvl-ipc.WINAC.SYSTEMSTATE.RED'**

**switchto\_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN\_REQUEST'**

**switchto\_operate = 'hvl-ipc.WINAC.SYSTEMSTATE.switchon'**

Writable boolean for switching to Red Operate (locket, HV on) state.

**switchto\_ready = 'hvl-ipc.WINAC.SYSTEMSTATE.RED\_REQUEST'**

Writable boolean for switching to Red Ready (locked, HV off) state.

**class SafetyStatus**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: [aenum.IntEnum](#)

Safety status values that are possible states returned from `hvl_ccb.dev.supercube.base.Supercube.get_status()`. These values correspond to the states of the Supercube’s safety circuit statemachine.

**Error = 6**

System is in error mode.

**GreenNotReady = 1**

System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

**GreenReady = 2**

System is safe and all safety elements are in place to be able to switch to *ready*.

**Initializing = 0**

System is initializing or booting.

**QuickStop = 5**

Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

**RedOperate = 4**

System is locked in red state and in *operate* mode, i.e. high voltage on.

**RedReady = 3**

System is locked in red state and *ready* to go to *operate* mode.

**class SupercubeOpcEndpoint**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl\_ccb.utils.enum.ValueEnum*

OPC Server Endpoint strings for the supercube variants.


**A = 'OPC.SimaticNET.S7'**

**B = 'OPC.SimaticNET.S7'**

**T13\_SOCKET\_PORTS = (1, 2, 3)**

Port numbers of SEV T13 power socket

**hvl\_ccb.dev.supercube2015.typ\_a**

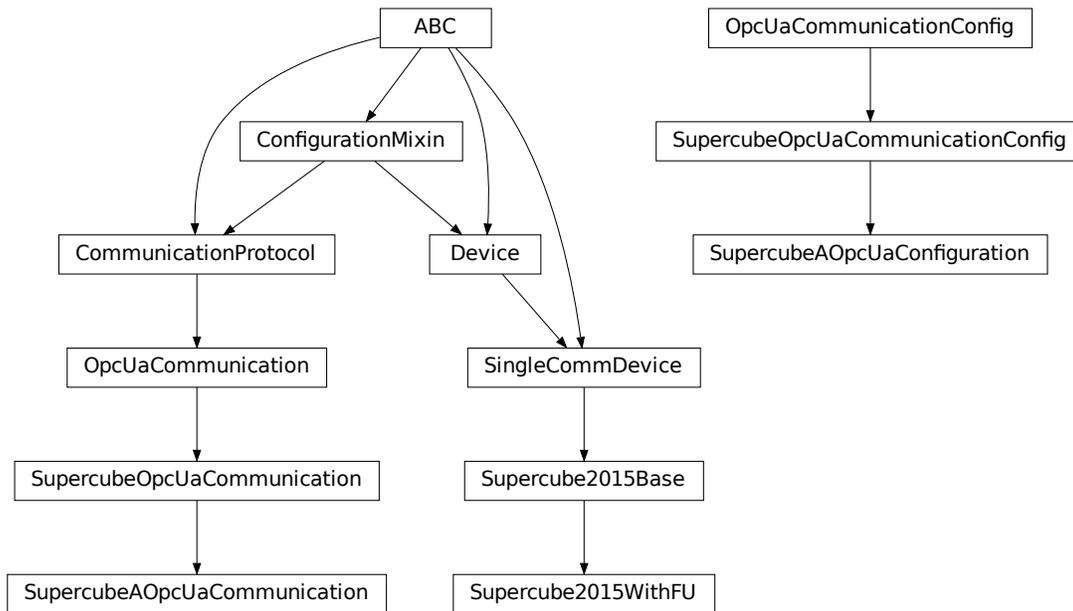

Supercube Typ A module.

**class Supercube2015WithFU**(*com*, *dev\_config=None*)

Bases: *hvl\_ccb.dev.supercube2015.base.Supercube2015Base*

Variant A of the Supercube with frequency converter.

**static default\_com\_cls**()

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**fso\_reset()** → *None*

Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.

**get\_frequency()** → float

Read the electrical frequency of the current Supercube setup.

**Returns** the frequency in Hz

**get\_fso\_active()** → bool

Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.

**Returns** state of the FSO functionality

**get\_max\_voltage()** → float

Reads the maximum voltage of the setup and returns in V.

**Returns** the maximum voltage of the setup in V.

**get\_power\_setup()** → *hvl\_ccb.dev.supercube2015.constants.PowerSetup*

Return the power setup selected in the Supercube's settings.

**Returns** the power setup

**get\_primary\_current()** → float

Read the current primary current at the output of the frequency converter ( before transformer).

**Returns** primary current in A

**get\_primary\_voltage()** → float

Read the current primary voltage at the output of the frequency converter ( before transformer).

**Returns** primary voltage in V

**get\_target\_voltage()** → float

Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to *set\_target\_voltage()*.

**Returns** the setpoint voltage in V.

**set\_slope(slope: float)** → *None*

Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.

**Parameters** **slope** – voltage slope in V/s (0..15'000)

**set\_target\_voltage(volt\_v: float)** → *None*

Set the output voltage to a defined value in V.

**Parameters** **volt\_v** – the desired voltage in V

**class SupercubeA0pcUaCommunication(config)**

Bases: *hvl\_ccb.dev.supercube2015.base.SupercubeOpcUaCommunication*

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class SupercubeA0pcUaConfiguration(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],
    endpoint_name: str = 'OPC.SimaticNET.S7', port: int = 4845,
    sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
    <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler
    object at 0x7f7a2c97d450>, update_parameter:
    asyncua.ua.uaprotoocol_auto.CreateSubscriptionParameters =
    CreateSubscriptionParameters(RequestedPublishingInterval=1000,
    RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22,
    MaxNotificationsPerPublish=10000, PublishingEnabled=True,
    Priority=0), wait_timeout_retry_sec: Union[int, float] = 1,
    max_timeout_retry_nr: int = 5)
```

Bases: `hvl_ccb.dev.supercube2015.base.SupercubeOpcUaCommunicationConfig`

**endpoint\_name:** `str = 'OPC.SimaticNET.S7'`

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


## Module contents

Supercube package with implementation for the old system version from 2015 based on Siemens WinAC soft-PLC on an industrial 32bit Windows computer.

`hvl_ccb.dev.technix`

Submodules

`hvl_ccb.dev.technix.base`

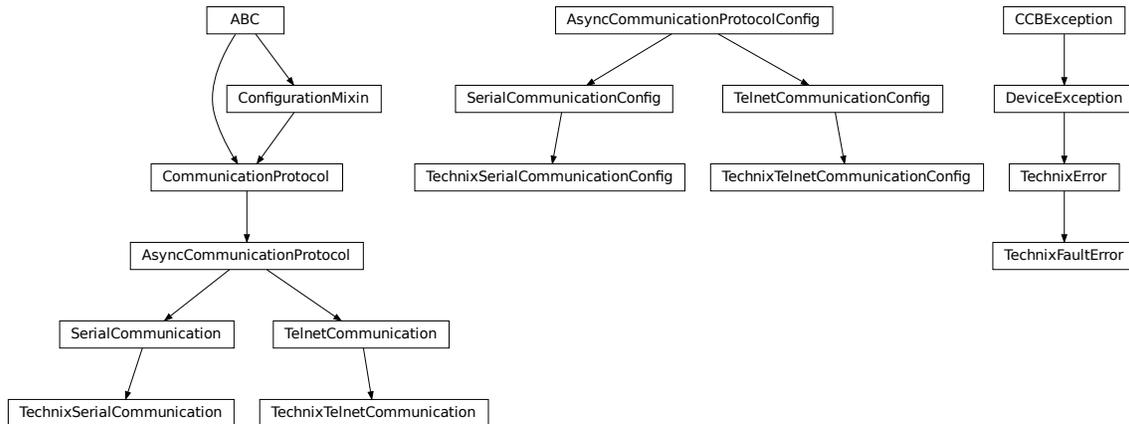

Communication and auxiliary classes for Technix

**exception TechnixError**

Bases: `hvl_ccb.dev.base.DeviceException`

Technix related errors.

**exception TechnixFaultError**

Bases: `hvl_ccb.dev.technix.base.TechnixError`

Raised when the fault flag was detected while the interlock is closed

**class TechnixSerialCommunication(configuration)**

Bases: `hvl_ccb.dev.technix.base._TechnixCommunication`, `hvl_ccb.comm.serial.SerialCommunication`

Serial communication for Technix

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class TechnixSerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                       encoding_error_handling: str = 'strict',
                                       wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                       default_n_attempts_read_text_nonempty: int = 10, port:
                                       Optional[str] = None, baudrate: int = 9600, parity: Union[str,
                                       hvl_ccb.comm.serial.SerialCommunicationParity] =
                                       SerialCommunicationParity.NONE, stopbits: Union[int, float,
                                       hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                       SerialCommunicationStopbits.ONE, bytesize: Union[int,
                                       hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                       SerialCommunicationBytesize.EIGHTBITS, timeout: Union[int,
                                       float] = 2)
```

Bases: `hvl_ccb.dev.technix.base._TechnixCommunicationConfig`, `hvl_ccb.comm.serial.SerialCommunicationConfig`

Configuration for the serial communication for Technix

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

```
class TechnixTelnetCommunication(configuration)
```

Bases: `hvl_ccb.comm.telnet.TelnetCommunication`, `hvl_ccb.dev.technix.base._TechnixCommunication`

Telnet communication for Technix

**static config\_cls**()

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class TechnixTelnetCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                     encoding_error_handling: str = 'strict',
                                     wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                     default_n_attempts_read_text_nonempty: int = 10, host:
                                     Optional[Union[str, ipaddress.IPv4Address,
                                     ipaddress.IPv6Address]] = None, port: int = 4660, timeout:
                                     Union[int, float] = 0.2)
```

Bases: `hvl_ccb.dev.technix.base._TechnixCommunicationConfig`, `hvl_ccb.comm.telnet.TelnetCommunicationConfig`

Configuration for the telnet communication for Technix

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 4660**


Port at which Technix is listening

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

## hvl\_ccb.dev.technix.device

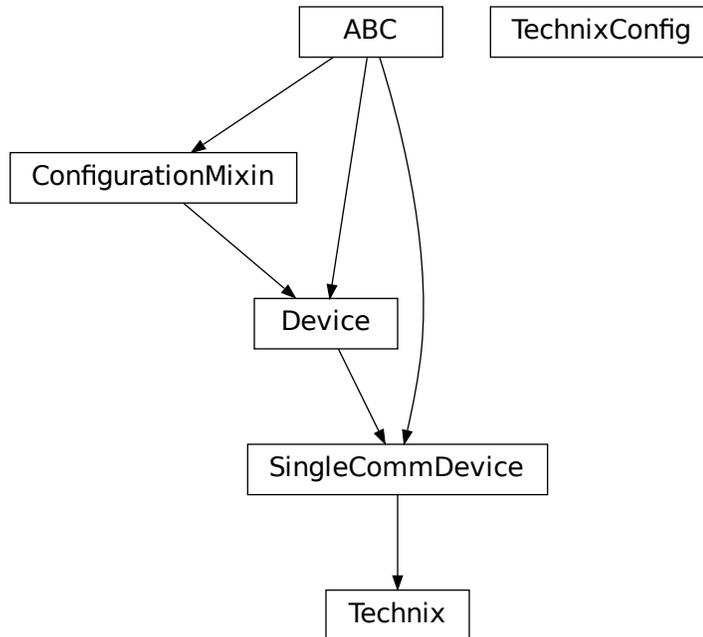

its corresponding configuration class

```
class Technix(com, dev_config)
```

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Device class to control capacitor chargers from Technix

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
property current: Optional[Union[int, float]]
```

Actual current of the output in A

```
default_com_cls() → Union[Type[hvl_ccb.dev.technix.base.TechnixSerialCommunication],  
Type[hvl_ccb.dev.technix.base.TechnixTelnetCommunication]]
```

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

```
property inhibit: Optional[bool]
```

Is the output of the voltage inhibited? The output stage can still be active.

```
property is_started: bool
```

Is the device started?

```
property max_current: Union[int, float]
```

Maximal output current of the hardware in A

The device class *Technix* and

**property max\_voltage: Union[int, float]**

Maximal output voltage of the hardware in V

**property open\_interlock: Optional[bool]**

Is the interlock open? (in safe mode)

**property output: Optional[bool]**

State of the high voltage output

**query\_status(\*, \_retry: bool = False)**

Query the status of the device.

**Returns** This function returns nothing

**property remote: Optional[bool]**

Is the device in remote control mode?

**start()**

Start the device and set it into the remote controllable mode. The high voltage is turn off, and the status poller is started.

**property status: Optional[hvl\_ccb.dev.technix.base.\_Status]**

The status of the device with the different states as sub-fields

**stop()**

Stop the device. The status poller is stopped and the high voltage output is turn off.

**property voltage: Optional[Union[int, float]]**

Actual voltage at the output in V

**property voltage\_regulation: Optional[bool]**

Status if the output is in voltage regulation mode (or current regulation)

**class TechnixConfig(communication\_channel:**

*Union[Type[hvl\_ccb.dev.technix.base.TechnixSerialCommunication],  
Type[hvl\_ccb.dev.technix.base.TechnixTelnetCommunication]], max\_voltage: Union[int,  
float], max\_current: Union[int, float], polling\_interval\_sec: Union[int, float] = 4,  
post\_stop\_pause\_sec: Union[int, float] = 1, register\_pulse\_time: Union[int, float] = 0.1,  
read\_output\_while\_polling: bool = False)*

Bases: object

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**communication\_channel:**

**Union[Type[hvl\_ccb.dev.technix.base.TechnixSerialCommunication],**

**Type[hvl\_ccb.dev.technix.base.TechnixTelnetCommunication]]**

communication channel between computer and Technix

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**max\_current: Union[int, float]**

Maximal Output current

**max\_voltage: Union[int, float]**

Maximal Output voltage

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_interval\_sec: Union[int, float] = 4**

Polling interval in s to maintain to watchdog of the device

**post\_stop\_pause\_sec: Union[int, float] = 1**

Time to wait after stopping the device

**read\_output\_while\_polling: bool = False**

Read output voltage and current within the polling event

**register\_pulse\_time: Union[int, float] = 0.1**

Time for pulsing a register

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

## Module contents

Device classes for “RS 232” and “Ethernet” interfaces, which are used to control power supplies from Technix. Manufacturer homepage: <https://www.technix-hv.com>

The regulated power supplies series and capacitor chargers series from Technix are series of low and high voltage direct current power supplies as well as capacitor chargers. The class *Technix* is tested with a CCR10KV-7,5KJ via an ethernet connection as well as a CCR15-P-2500-OP via a serial connection. Check the code carefully before using it with other devices or device series


**This Python package may support the following interfaces from Technix:**

- [Remote Interface RS232](#)
- [Ethernet Remote Interface](#)
- [Optic Fiber Remote Interface](#)

`hvl_ccb.dev.tiepie`

Submodules

`hvl_ccb.dev.tiepie.base`

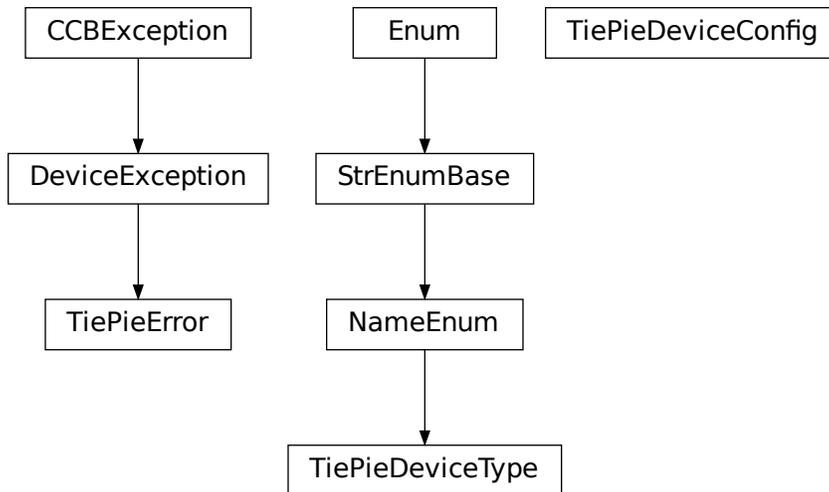

```

class TiePieDeviceConfig(serial_number: int, require_block_measurement_support: bool = True,
                        n_max_try_get_device: int = 10, wait_sec_retry_get_device: Union[int, float] =
                        1.0, is_data_ready_polling_interval_sec: Union[int, float] = 0.01)
  
```

Bases: object

Configuration dataclass for TiePie

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**is\_data\_ready\_polling\_interval\_sec: Union[int, float] = 0.01**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**n\_max\_try\_get\_device:** `int = 10`

**classmethod optional\_defaults()** `→ Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**require\_block\_measurement\_support:** `bool = True`

**classmethod required\_keys()** `→ Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**serial\_number:** `int`

**wait\_sec\_retry\_get\_device:** `Union[int, float] = 1.0`

**class TiePieDeviceType** (*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `hvl_ccb.utils.enum.NameEnum`

TiePie device type.

**GENERATOR = 2**

**I2C = 4**

**OSCILLOSCOPE = 1**

**exception TiePieError**

Bases: `hvl_ccb.dev.base.DeviceException`

Error of the class TiePie

**get\_device\_by\_serial\_number** (*serial\_number: int, device\_type: Union[str, Tuple[int, hvl\_ccb.dev.tiepie.base.LtpDeviceReturnType]], n\_max\_try\_get\_device: int = 10, wait\_sec\_retry\_get\_device: float = 1.0*) `→ hvl_ccb.dev.tiepie.base.LtpDeviceReturnType`

Open and return handle of TiePie device with a given serial number

#### Parameters

- **serial\_number** – int serial number of the device
- **device\_type** – a `TiePieDeviceType` instance containing device identifier (int number) and its corresponding class, both from `libtiepie`, or a string name of such instance
- **n\_max\_try\_get\_device** – maximal number of device list updates (int number)
- **wait\_sec\_retry\_get\_device** – waiting time in seconds between retries (int number)

**Returns** Instance of a `libtiepie` device class according to the specified `device_type`

#### Raises

- **TiePieError** – when there is no device with given serial number
- **ValueError** – when `device_type` is not an instance of `TiePieDeviceType`

**wrap\_libtiepie\_exception** (*func: Callable*) `→ Callable`


Decorator wrapper for `libtiepie` methods that use `libtiepie.library.check_last_status_raise_on_error()` calls.

**Parameters** **func** – Function or method to be wrapped

**Raises** *TiePieError* – instead of *LibTiePieException* or one of its subtypes.

**Returns** whatever *func* returns

## hvl\_ccb.dev.tiepie.channel

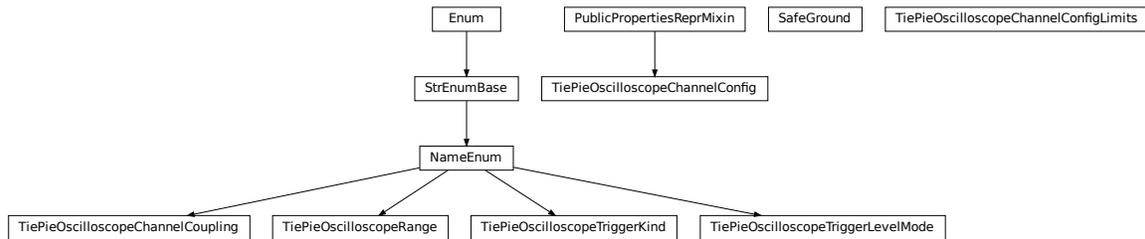

### class SafeGround

Bases: object

Class that dynamically adds the `safe_ground_enabled` attribute getter/setter if the bound oscilloscope has the safe ground option.

### class TiePieOscilloscopeChannelConfig(*ch\_number: int, channel: libtiepie.oscilloscopechannel.OscilloscopeChannel*)

Bases: *hvl\_ccb.dev.tiepie.utils.PublicPropertiesReprMixin*

Oscilloscope's channel configuration, with cleaning of values in properties setters as well as setting and reading them on and from the device's channel.

**static clean\_coupling**(*coupling: Union[str, hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling]*) → *hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling*

**static clean\_enabled**(*enabled: bool*) → bool

**clean\_input\_range**(*input\_range: Union[float, hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeRange]*) → *hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeRange*

**clean\_probe\_offset**(*probe\_offset: float*) → float

**static clean\_trigger\_enabled**(*trigger\_enabled*)

**clean\_trigger\_hysteresis**(*trigger\_hysteresis: float*) → float

**static clean\_trigger\_kind**(*trigger\_kind: Union[str, hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind]*) → *hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind*

**clean\_trigger\_level**(*trigger\_level: Union[int, float]*) → float

**static clean\_trigger\_level\_mode**(*level\_mode: Union[str, hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode]*) → *hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode*

property `coupling`: *hvl\_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling*

property `enabled`: bool

**property has\_safe\_ground: bool**

Check whether bound oscilloscope device has “safe ground” option

**Returns** bool: 1=safe ground available

**property input\_range:** [hvl\\_ccb.dev.tiepie.channel.TiePieOscilloscopeRange](#)

**property probe\_offset:** float

**property trigger\_enabled:** bool

**property trigger\_hysteresis:** float

**property trigger\_kind:** [hvl\\_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind](#)

**property trigger\_level:** float

**property trigger\_level\_mode:**

[hvl\\_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode](#)

**class TiePieOscilloscopeChannelConfigLimits**(*osc\_channel:*  
*libtiepie.oscilloscopechannel.OscilloscopeChannel*)

Bases: object

Default limits for oscilloscope channel parameters.

**class TiePieOscilloscopeChannelCoupling**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*,  
*start=1*, *boundary=None*)

Bases: [hvl\\_ccb.utils.enum.NameEnum](#)

An enumeration.

**ACA = 8**

**ACV = 2**

**DCA = 4**

**DCV = 1**

**class TiePieOscilloscopeRange**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*,  
*boundary=None*)

Bases: [hvl\\_ccb.utils.enum.NameEnum](#)

An enumeration.

**EIGHTY\_VOLT = 80**

**EIGHT\_HUNDRED\_MILLI\_VOLT = 0.8**

**EIGHT\_VOLT = 8**

**FORTY\_VOLT = 40**

**FOUR\_HUNDRED\_MILLI\_VOLT = 0.4**

**FOUR\_VOLT = 4**

**TWENTY\_VOLT = 20**

**TWO\_HUNDRED\_MILLI\_VOLT = 0.2**

**TWO\_VOLT = 2**

**static suitable\_range**(*value*)

**class TiePieOscilloscopeTriggerKind**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl\_ccb.utils.enum.NameEnum*

An enumeration.

**ANY** = 16

**FALLING** = 2

**RISING** = 1

**RISING\_OR\_FALLING** = 16

**class TiePieOscilloscopeTriggerLevelMode**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl\_ccb.utils.enum.NameEnum*


An enumeration.

**ABSOLUTE** = 2

**RELATIVE** = 1

**UNKNOWN** = 0

**hvl\_ccb.dev.tiepie.device**

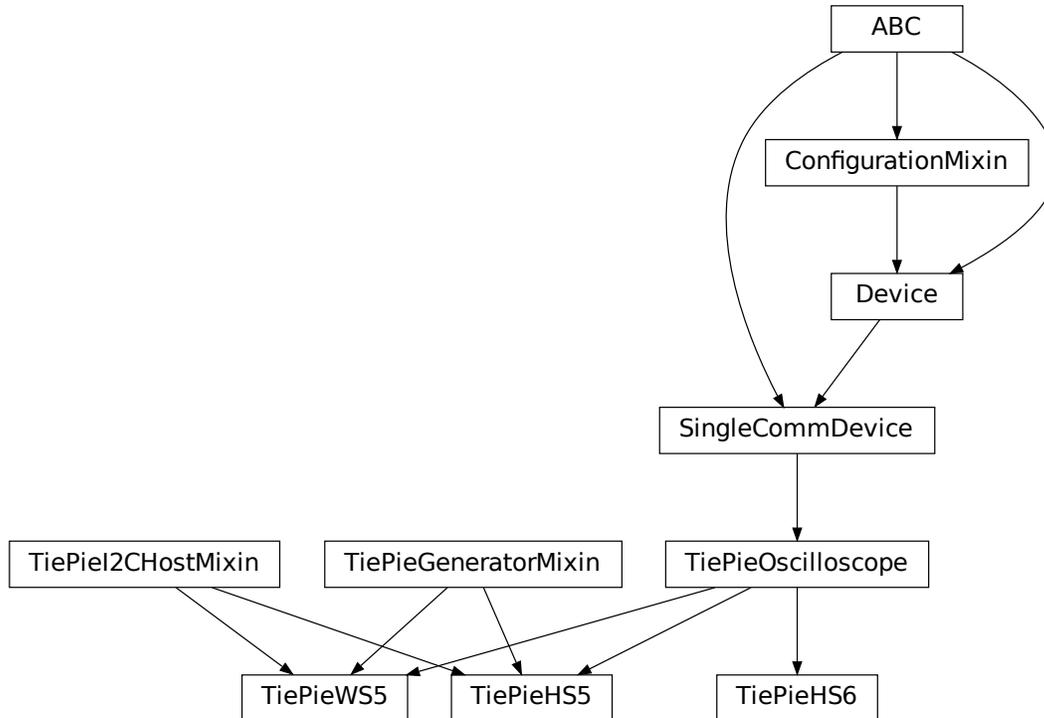

TiePie devices.

```
class TiePieHS5(com, dev_config)
```

```
Bases: hvl_ccb.dev.tiepie.i2c.TiePieI2CHostMixin, hvl_ccb.dev.tiepie.generator.TiePieGeneratorMixin, hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope
```

TiePie HS5 device.

```
class TiePieHS6(com, dev_config)
```

```
Bases: hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope
```


TiePie HS6 DIFF device.

```
class TiePieWS5(com, dev_config)
```

```
Bases: hvl_ccb.dev.tiepie.i2c.TiePieI2CHostMixin, hvl_ccb.dev.tiepie.generator.TiePieGeneratorMixin, hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope
```

TiePie WS5 device.

### **hvl\_ccb.dev.tiepie.generator**

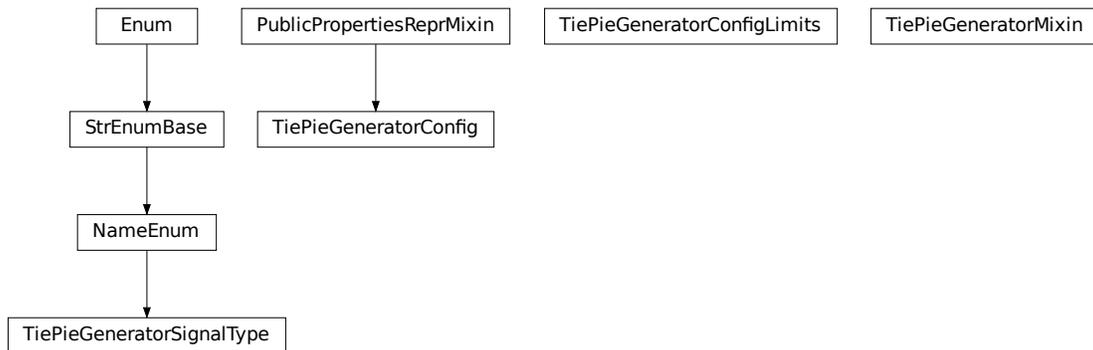

```
class TiePieGeneratorConfig(dev_gen: libtiepie.generator.Generator)
```

```
Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin
```

Generator's configuration with cleaning of values in properties setters.

```
property amplitude: float
```

```
clean_amplitude(amplitude: float) → float
```

```
static clean_enabled(enabled: bool) → bool
```

```
clean_frequency(frequency: float) → float
```

```
clean_offset(offset: float) → float
```

```
static clean_signal_type(signal_type: Union[int,  
hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType]) →  
hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType
```

```
clean_waveform(waveform: numpy.ndarray[Any, numpy.dtype[numpy.typing._generic_alias.ScalarType]])  
→ numpy.ndarray[Any, numpy.dtype[numpy.typing._generic_alias.ScalarType]]
```

```
property enabled: bool
```

property frequency: float

property offset: float

property signal\_type: *hvl\_ccb.dev.tiepie.generator.TiePieGeneratorSignalType*

property waveform: Optional[numpy.ndarray[Any,  
numpy.dtype[numpy.typing.\_generic\_alias.ScalarType]]]

**class TiePieGeneratorConfigLimits**(*dev\_gen: libtiepie.generator.Generator*)

Bases: object

Default limits for generator parameters.

**class TiePieGeneratorMixin**(*com, dev\_config*)

Bases: object

TiePie Generator sub-device.

A wrapper for the *libtiepie.generator.Generator* class. To be mixed in with *TiePieOscilloscope* base class.

**config\_gen:** Optional[*hvl\_ccb.dev.tiepie.generator.TiePieGeneratorConfig*]

Generator's dynamical configuration.

**generator\_start()**

Start signal generation.

**generator\_stop()**

Stop signal generation.

**start()** → *None*

Start the Generator.

**stop()** → *None*

Stop the generator.

**class TiePieGeneratorSignalType**(*value=<no\_arg>, names=None, module=None, type=None, start=1,  
boundary=None*)

Bases: *hvl\_ccb.utils.enum.NameEnum*

An enumeration.

**ARBITRARY = 32**

**DC = 8**

**NOISE = 16**

**PULSE = 64**


**SINE = 1**

**SQUARE = 4**

**TRIANGLE = 2**

**UNKNOWN = 0**

## hvl\_ccb.dev.tiepie.i2c

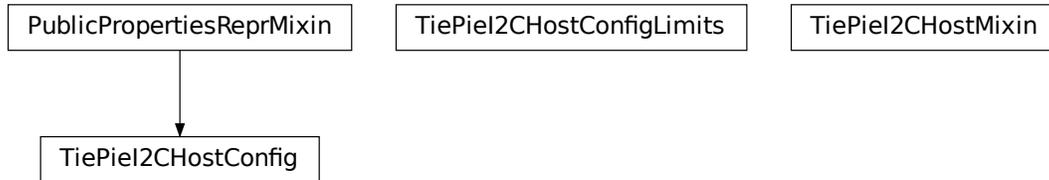

**class TiePieI2CHostConfig**(*dev\_i2c: libtiepie.i2chost.I2CHost*)

Bases: *hvl\_ccb.dev.tiepie.utils.PublicPropertiesReprMixin*

I2C Host's configuration with cleaning of values in properties setters.

**class TiePieI2CHostConfigLimits**(*dev\_i2c: libtiepie.i2chost.I2CHost*)

Bases: object

Default limits for I2C host parameters.

**class TiePieI2CHostMixin**(*com, dev\_config*)

Bases: object

TiePie I2CHost sub-device.

A wrapper for the *libtiepie.i2chost.I2CHost* class. To be mixed in with *TiePieOscilloscope* base class.

**config\_i2c: Optional**[*hvl\_ccb.dev.tiepie.i2c.TiePieI2CHostConfig*]

I2C host's dynamical configuration.


**start()** → *None*

Start the I2C Host.

**stop()** → *None*

Stop the I2C host.

## hvl\_ccb.dev.tiepie.oscilloscope

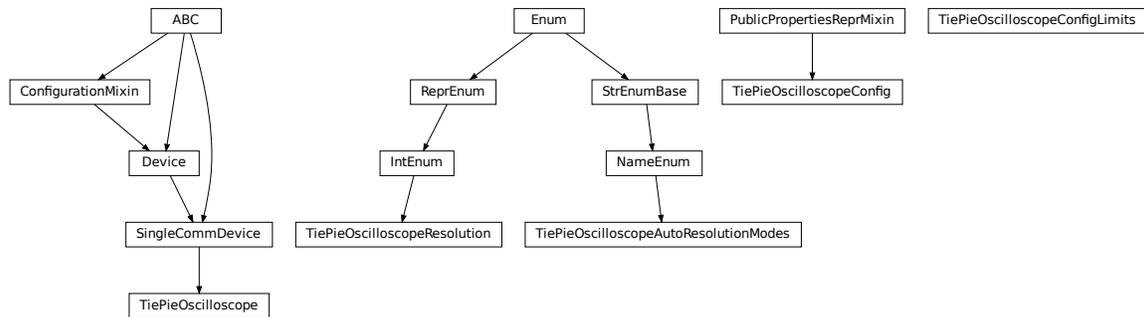

**class** `TiePieOscilloscope`(*com*, *dev\_config*)

Bases: `hvl_ccb.dev.base.SingleCommDevice`

TiePie oscilloscope.

A wrapper for TiePie oscilloscopes, based on the class `libtiepie.oscilloscope.Oscilloscope` with simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and its channels, including extra validation and typing hints support for configurations.

Note that, in contrast to `libtiepie` library, since all physical TiePie devices include an oscilloscope, this is the base class for all physical TiePie devices. The additional TiePie sub-devices: “Generator” and “I2CHost”, are mixed-in to this base class in subclasses.

The channels use  $1..N$  numbering (not  $0..N-1$ ), as in, e.g., the Multi Channel software.

**property** `channels_enabled`: `Generator[int, None, None]`

Yield numbers of enabled channels.

**Returns** Numbers of enabled channels

**collect\_measurement\_data**(*timeout*: `Optional[Union[int, float]] = 0`) → `Optional[numpy.ndarray[Any, numpy.dtype[numpy.typing._generic_alias.ScalarType]]]`

Try to collect the data from TiePie; return `None` if data is not ready.

**Parameters** `timeout` – The timeout to wait until data is available. This option makes this function blocking the code. `timeout = None` blocks the code infinitely till data will be available. Per default, the `timeout` is set to `0`: The function will not block.

**Returns** Measurement data of only enabled channels and time vector in a `2D-numpy.ndarray` with float sample data; or `None` if there is no data available.

**static** `config_cls`() → `Type[hvl_ccb.dev.tiepie.base.TiePieDeviceConfig]`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**config\_osc**: `Optional[hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeConfig]`

Oscilloscope’s dynamical configuration.

**config\_osc\_channel\_dict**: `Dict[int, hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelConfig]`

Channel configuration. A `dict` mapping actual channel number, numbered  $1..N$ , to channel configuration. The channel info is dynamically read from the device only on the first `start()`; beforehand the `dict` is empty.

**static default\_com\_cls()** → Type[hvl\_ccb.comm.base.NullCommunicationProtocol]

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**force\_trigger()** → None

Forces the TiePie to trigger with a software sided trigger event.

**Return** None

**Raises** *TiePieError* – when device is not started or status of underlying device gives an error

**is\_measurement\_data\_ready()** → bool

Reports if TiePie has data which is ready to collect

**Returns** if the data is ready to collect.

**Raises** *TiePieError* – when device is not started or status of underlying device gives an error

**is\_triggered()** → bool

Reports if TiePie has triggered. Maybe data is not yet available. One can check with the function *is\_measurement\_data\_ready()*.

**Returns** if a trigger event occurred

**static list\_devices()** → libtiepie.devicelist.DeviceList

List available TiePie devices.

**Returns** libtiepie up to date list of devices

**property n\_channels**

Number of channels in the oscilloscope.

**Returns** Number of channels.

**start()** → None

Start the oscilloscope.

**start\_measurement()** → None

Start a measurement using set configuration.

**Raises** *TiePieError* – when device is not started or status of underlying device gives an error

**stop()** → None

Stop the oscilloscope.

**class TiePieOscilloscopeAutoResolutionModes**(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: *hvl\_ccb.utils.enum.NameEnum*

An enumeration.

**ALL** = 4

**DISABLED** = 1

**NATIVEONLY** = 2

**UNKNOWN** = 0

**class TiePieOscilloscopeConfig**(dev\_osc: libtiepie.oscilloscope.Oscilloscope)

Bases: *hvl\_ccb.dev.tiepie.utils.PublicPropertiesReprMixin*

Oscilloscope's configuration with cleaning of values in properties setters.

**property auto\_resolution\_mode:**

*hvl\_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes*

```
static clean_auto_resolution_mode(auto_resolution_mode: Union[int,
                                hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes])
    →
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes

clean_pre_sample_ratio(pre_sample_ratio: float) → float
clean_record_length(record_length: Union[int, float]) → int
static clean_resolution(resolution: Union[int,
                                           hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution]) →
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution

clean_sample_frequency(sample_frequency: float) → float
clean_trigger_timeout(trigger_timeout: Optional[Union[int, float]]) → float
property pre_sample_ratio: float
property record_length: int
property resolution: hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution
property sample_frequency: float
property trigger_timeout: Optional[float]
```

```
class TiePieOscilloscopeConfigLimits(dev_osc: libtiepie.oscilloscope.Oscilloscope)
```

Bases: object

Default limits for oscilloscope parameters.

```
class TiePieOscilloscopeResolution(value=<no_arg>, names=None, module=None, type=None, start=1,
                                   boundary=None)
```

Bases: aenum.IntEnum

An enumeration.

**EIGHT\_BIT = 8**

**FOURTEEN\_BIT = 14**

**SIXTEEN\_BIT = 16**

**TWELVE\_BIT = 12**

## [hvl\\_ccb.dev.tiepie.utils](#)

PublicPropertiesReprMixin

```
class PublicPropertiesReprMixin
```

Bases: object

General purpose utility mixin that overwrites object representation to a one analogous to *dataclass* instances, but using public properties and their values instead of *fields*.

## Module contents

This module is a wrapper around LibTiePie SDK devices; see <https://www.tiepie.com/en/libtiepie-sdk> .

The device classes adds simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and oscilloscope's channels. This includes extra validation and typing hints support.

## Extra installation

LibTiePie SDK library is available only on Windows and on Linux.

To use this LibTiePie SDK devices wrapper:

1. install the `hvl_ccb` package with a `tiepie` extra feature:

```
$ pip install "hvl_ccb[tiepie]"
```

this will install the Python bindings for the library.

2. install the library

- on Linux: follow instructions in <https://www.tiepie.com/en/libtiepie-sdk/linux> ;
- on Windows: the additional DLL is included in Python bindings package.

## Troubleshooting

On a Windows system, if you encounter an `OSError` like this:

```
...
    self._handle = _dlopen(self._name, mode)
OSError: [WinError 126] The specified module could not be found
```


most likely the `python-libtiepie` package was installed in your `site-packages/` directory as a `python-libtiepie-*.egg` file via `python setup.py install` or `python setup.py develop` command. In such case uninstall the library and re-install it using `pip`:

```
$ pip uninstall python-libtiepie
$ pip install python-libtiepie
```

This should create `libtiepie/` folder. Alternatively, manually move the folder `libtiepie/` from inside of the `.egg` archive file to the containing it `site-packages/` directory (PyCharm's Project tool window supports reading and extracting from `.egg` archives).

## Submodules

## hvl\_ccb.dev.base

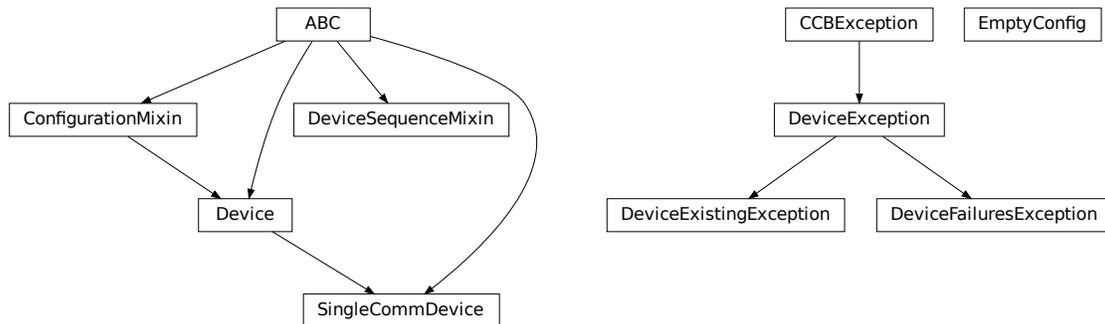

Module with base classes for devices.

**class** `Device`(*dev\_config=None*)

Bases: `hvl_ccb.configuration.ConfigurationMixin`, `abc.ABC`

Base class for devices. Implement this class for a concrete device, such as measurement equipment or voltage sources.

Specifies the methods to implement for a device.

**static** `config_cls()`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**abstract** `start()` → *None*

Start or restart this Device. To be implemented in the subclass.

**abstract** `stop()` → *None*

Stop this Device. To be implemented in the subclass.

**exception** `DeviceException`

Bases: `hvl_ccb.exception.CCBException`

**exception** `DeviceExistingException`

Bases: `hvl_ccb.dev.base.DeviceException`

Exception to indicate that a device with that name already exists.

**exception** `DeviceFailuresException`(*failures: Dict[str, Exception], \*args*)

Bases: `hvl_ccb.dev.base.DeviceException`

Exception to indicate that one or several devices failed.

**failures:** `Dict[str, Exception]`

A dictionary of named devices failures (exceptions).

**class** `DeviceSequenceMixin`(*devices: Dict[str, hvl\_ccb.dev.base.Device]*)

Bases: `abc.ABC`

Mixin that can be used on a device or other classes to provide facilities for handling multiple devices in a sequence.

**add\_device**(*name: str, device: hvl\_ccb.dev.base.Device*) → *None*

Add a new device to the device sequence.

**Parameters**

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

**Raises** *DeviceExistingException* –

**devices\_failed\_start:** Dict[str, *hvl\_ccb.dev.base.Device*]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices\_failed\_stop*.

**devices\_failed\_stop:** Dict[str, *hvl\_ccb.dev.base.Device*]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices\_failed\_start*.

**get\_device**(*name: str*) → *hvl\_ccb.dev.base.Device*

Get a device by name.

**Parameters** **name** – is the name of the device.

**Returns** the device object from this sequence.

**get\_devices**() → List[Tuple[str, *hvl\_ccb.dev.base.Device*]]

Get list of name, device pairs according to current sequence.

**Returns** A list of tuples with name and device each.

**remove\_device**(*name: str*) → *hvl\_ccb.dev.base.Device*

Remove a device from this sequence and return the device object.

**Parameters** **name** – is the name of the device.

**Returns** device object or *None* if such device was not in the sequence.

**Raises** **ValueError** – when device with given name was not found

**start**() → *None*

Start all devices in this sequence in their added order.

**Raises** *DeviceFailuresException* – if one or several devices failed to start

**stop**() → *None*

Stop all devices in this sequence in their reverse order.

**Raises** *DeviceFailuresException* – if one or several devices failed to stop

**class** **EmptyConfig**

Bases: object

Empty configuration dataclass that is the default configuration for a Device.

**clean\_values**()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**class SingleCommDevice**(*com, dev\_config=None*)

Bases: *hvl\_ccb.dev.base.Device*, *abc.ABC*

Base class for devices with a single communication protocol.

**property com**

Get the communication protocol of this device.

**Returns** an instance of CommunicationProtocol subtype

**abstract static default\_com\_cls()** → Type[*hvl\_ccb.comm.base.CommunicationProtocol*]

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device


**start()** → *None*

Open the associated communication protocol.

**stop()** → *None*

Close the associated communication protocol.

## hvl\_ccb.dev.crylas

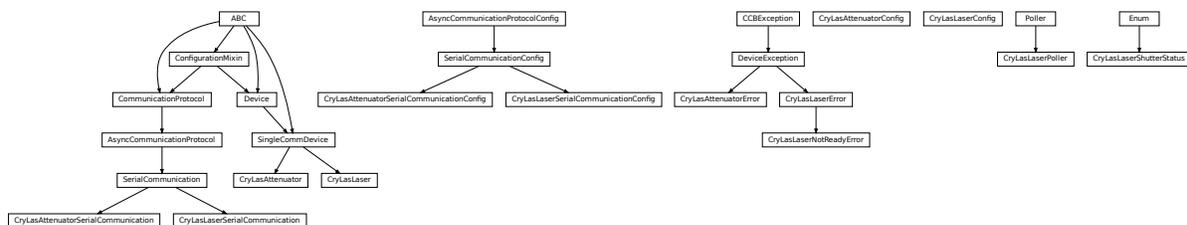

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: [https://www.crylas.de/products/pulsed\\_laser.html](https://www.crylas.de/products/pulsed_laser.html)

**class CryLasAttenuator**(*com, dev\_config=None*)

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Device class for the CryLas laser attenuator.

**property attenuation:** `Union[int, float]`

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**set\_attenuation**(*percent: Union[int, float]*) → *None*

Set the percentage of attenuated light (inverse of `set_transmission`). :param *percent*: percentage of attenuation, number between 0 and 100 :raises `ValueError`: if param *percent* not between 0 and 100 :raises `SerialCommunicationIOError`: when communication port is not opened :raises `CryLasAttenuatorError`: if the device does not confirm success

**set\_init\_attenuation()**

Sets the attenuation to its configured initial/default value

**Raises** `SerialCommunicationIOError` – when communication port is not opened

**set\_transmission**(*percent: Union[int, float]*) → *None*

Set the percentage of transmitted light (inverse of `set_attenuation`). :param *percent*: percentage of transmitted light :raises `ValueError`: if param *percent* not between 0 and 100 :raises `SerialCommunicationIOError`: when communication port is not opened :raises `CryLasAttenuatorError`: if the device does not confirm success

**start()** → *None*

Open the com, apply the config value 'init\_attenuation'

**Raises** `SerialCommunicationIOError` – when communication port cannot be opened

**property transmission:** `Union[int, float]`

**class CryLasAttenuatorConfig**(*init\_attenuation: Union[int, float] = 0, response\_sleep\_time: Union[int, float] = 1*)

Bases: `object`

Device configuration dataclass for CryLas attenuator.

**clean\_values()**

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**init\_attenuation:** `Union[int, float] = 0`

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**response\_sleep\_time: Union[int, float] = 1**

**exception CryLasAttenuatorError**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

General error with the CryLas Attenuator.

**class CryLasAttenuatorSerialCommunication(configuration)**

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for the CryLas attenuator. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class CryLasAttenuatorSerialCommunicationConfig(terminator: bytes = b'', encoding: str = 'utf-8',
                                                encoding_error_handling: str = 'strict',
                                                wait_sec_read_text_nonempty: Union[int, float] =
                                                0.5, default_n_attempts_read_text_nonempty: int =
                                                10, port: Union[str, NoneType] = None, baudrate: int
                                                = 9600, parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize]
                                                = <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 3)
```

Bases: [hvl\\_ccb.comm.serial.SerialCommunicationConfig](#)

**baudrate: int = 9600**

Baudrate for CryLas attenuator is 9600 baud

**bytesize: Union[int, [hvl\\_ccb.comm.serial.SerialCommunicationBytesize](#)] = 8**

One byte is eight bits long

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** Union[str, [hvl\\_ccb.comm.serial.SerialCommunicationParity](#)] = 'N'  
CryLas attenuator does not use parity

**classmethod required\_keys()** → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** Union[int, [hvl\\_ccb.comm.serial.SerialCommunicationStopbits](#)] = 1  
CryLas attenuator uses one stop bit

**terminator:** bytes = b''  
No terminator

**timeout:** Union[int, float] = 3  
use 3 seconds timeout as default

**class CryLasLaser**(*com*, *dev\_config=None*)  
Bases: [hvl\\_ccb.dev.base.SingleCommDevice](#)

CryLas laser controller device class.

**class AnswersShutter**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [aenum.Enum](#)

Standard answers of the CryLas laser controller to 'Shutter' command passed via *com*.

**CLOSED** = 'Shutter inaktiv'

**OPENED** = 'Shutter aktiv'

**class AnswersStatus**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [aenum.Enum](#)

Standard answers of the CryLas laser controller to 'STATUS' command passed via *com*.

**ACTIVE** = 'STATUS: Laser active'

**HEAD** = 'STATUS: Head ok'

**INACTIVE** = 'STATUS: Laser inactive'

**READY** = 'STATUS: System ready'

```
TEC1 = 'STATUS: TEC1 Regulation ok'
TEC2 = 'STATUS: TEC2 Regulation ok'
```

**class LaserStatus**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)  
Bases: aenum.Enum  
Status of the CryLas laser

```
READY_ACTIVE = 2
READY_INACTIVE = 1
UNREADY_INACTIVE = 0
property is_inactive
property is_ready
```

**class RepetitionRates**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*,  
*boundary=None*)  
Bases: aenum.IntEnum  
Repetition rates for the internal software trigger in Hz

```
HARDWARE = 0
SOFTWARE_INTERNAL_SIXTY = 60
SOFTWARE_INTERNAL_TEN = 10
SOFTWARE_INTERNAL_TWENTY = 20
```

**ShutterStatus**  
alias of *hvl\_ccb.dev.crylas.CryLasLaserShutterStatus*

**close\_shutter()** → *None*  
Close the laser shutter.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**static config\_cls()**  
Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**  
Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**get\_pulse\_energy\_and\_rate()** → Tuple[int, int]  
Use the debug mode, return the measured pulse energy and rate.

**Returns** (energy in micro joule, rate in Hz)

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not answer the query

**laser\_off()** → *None*  
Turn the laser off.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**laser\_on()** → *None*

Turn the laser on.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserNotReadyError* – if the laser is not ready to be turned on
- *CryLasLaserError* – if success is not confirmed by the device

**open\_shutter()** → *None*

Open the laser shutter.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**set\_init\_shutter\_status()** → *None*

Open or close the shutter, to match the configured shutter\_status.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**set\_pulse\_energy(energy: int)** → *None*

Sets the energy of pulses (works only with external hardware trigger). Proceed with small energy steps, or the regulation may fail.

**Parameters energy** – energy in micro joule

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not confirm success

**set\_repetition\_rate(rate: Union[int, hvl\_ccb.dev.crylas.CryLasLaser.RepetitionRates])** → *None*

Sets the repetition rate of the internal software trigger.

**Parameters rate** – frequency (Hz) as an integer

**Raises**

- **ValueError** – if rate is not an accepted value in RepetitionRates Enum
- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**start()** → *None*

Opens the communication protocol and configures the device.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened

**stop()** → *None*

Stops the device and closes the communication protocol.

**Raises**

- ***SerialCommunicationIOError*** – if com port is closed unexpectedly
- ***CryLasLaserError*** – if `laser_off()` or `close_shutter()` fail

**property** `target_pulse_energy`

**update\_laser\_status()** → *None*

Update the laser status to *LaserStatus.NOT\_READY* or *LaserStatus.INACTIVE* or *LaserStatus.ACTIVE*.

Note: laser never explicitly says that it is not ready (*LaserStatus.NOT\_READY*) in response to ‘STATUS’ command. It only says that it is ready (heated-up and implicitly inactive/off) or active (on). If it’s not either of these then the answer is *Answers.HEAD*. Moreover, the only time the laser explicitly says that its status is inactive (*Answers.INACTIVE*) is after issuing a ‘LASER OFF’ command.

**Raises** ***SerialCommunicationIOError*** – when communication port is not opened

**update\_repetition\_rate()** → *None*

Query the laser repetition rate.

**Raises**

- ***SerialCommunicationIOError*** – when communication port is not opened
- ***CryLasLaserError*** – if success is not confirmed by the device

**update\_shutter\_status()** → *None*

Update the shutter status (OPENED or CLOSED)

**Raises**

- ***SerialCommunicationIOError*** – when communication port is not opened
- ***CryLasLaserError*** – if success is not confirmed by the device

**update\_target\_pulse\_energy()** → *None*

Query the laser pulse energy.

**Raises**

- ***SerialCommunicationIOError*** – when communication port is not opened
- ***CryLasLaserError*** – if success is not confirmed by the device

**wait\_until\_ready()** → *None*

Block execution until the laser is ready

**Raises** ***CryLasLaserError*** – if the polling thread stops before the laser is ready

```
class CryLasLaserConfig(calibration_factor: Union[int, float] = 4.35, polling_period: Union[int, float] = 12,  
                       polling_timeout: Union[int, float] = 300, auto_laser_on: bool = True,  
                       init_shutter_status: Union[int, hvl_ccb.dev.crylas.CryLasLaserShutterStatus] =  
                       CryLasLaserShutterStatus.CLOSED)
```

Bases: object

Device configuration dataclass for the CryLas laser controller.

**ShutterStatus**

alias of *hvl\_ccb.dev.crylas.CryLasLaserShutterStatus*

**auto\_laser\_on:** bool = True

**calibration\_factor:** Union[int, float] = 4.35

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**init\_shutter\_status:** Union[int, [hvl\\_ccb.dev.crylas.CryLasLaserShutterStatus](#)] = 0

**is\_configdataclass** = True

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_period:** Union[int, float] = 12

**polling\_timeout:** Union[int, float] = 300

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception CryLasLaserError**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

General error with the CryLas Laser.

**exception CryLasLaserNotReadyError**

Bases: [hvl\\_ccb.dev.crylas.CryLasLaserError](#)

Error when trying to turn on the CryLas Laser before it is ready.

**class CryLasLaserPoller**(*spoll\_handler*: Callable, *check\_handler*: Callable, *check\_laser\_status\_handler*: Callable, *polling\_delay\_sec*: Union[int, float] = 0, *polling\_interval\_sec*: Union[int, float] = 1, *polling\_timeout\_sec*: Optional[Union[int, float]] = None)

Bases: [hvl\\_ccb.dev.utils.Poller](#)

Poller class for polling the laser status until the laser is ready.

**Raises**

- **CryLasLaserError** – if the timeout is reached before the laser is ready
- **SerialCommunicationIOError** – when communication port is closed.

**class CryLasLaserSerialCommunication**(*configuration*)

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for the CryLas laser controller. Already predefines device-specific protocol parameters in config.

**READ\_TEXT\_SKIP\_PREFIXES** = ('>', 'MODE:')

Prefixes of lines that are skipped when read from the serial port.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query**(*cmd: str, prefix: str, post\_cmd: Optional[str] = None*) → str

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns the line in question.

**Parameters**

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer
- **post\_cmd** – optional additional command to send after the query

**Returns** line in question as a string

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**query\_all**(*cmd: str, prefix: str*)

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns a list of successive lines starting with prefix.

**Parameters**

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer

**Returns** line in question as a string

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**read**() → str

Read first line of text from the serial port that does not start with any of *self.READ\_TEXT\_SKIP\_PREFIXES*.

**Returns** String read from the serial port; '' if there was nothing to read.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

```
class CryLasLaserSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',  
                                           encoding_error_handling: str = 'strict',  
                                           wait_sec_read_text_nonempty: Union[int, float] = 0.5,  
                                           default_n_attempts_read_text_nonempty: int = 10, port:  
                                           Union[str, NoneType] = None, baudrate: int = 19200,  
                                           parity: Union[str,  
                                           hvl_ccb.comm.serial.SerialCommunicationParity] =  
                                           <SerialCommunicationParity.NONE: 'N'>, stopbits:  
                                           Union[int,  
                                           hvl_ccb.comm.serial.SerialCommunicationStopbits] =  
                                           <SerialCommunicationStopbits.ONE: 1>, bytesize:  
                                           Union[int,  
                                           hvl_ccb.comm.serial.SerialCommunicationBytesize] =  
                                           <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:  
                                           Union[int, float] = 10)
```

Bases: *hvl\_ccb.comm.serial.SerialCommunicationConfig*

**baudrate: int = 19200**

Baudrate for CryLas laser is 19200 baud

**bytesize: Union[int, *hvl\_ccb.comm.serial.SerialCommunicationBytesize*] = 8**

One byte is eight bits long

**force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity: Union[str, *hvl\_ccb.comm.serial.SerialCommunicationParity*] = 'N'**

CryLas laser does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits: Union[int, *hvl\_ccb.comm.serial.SerialCommunicationStopbits*] = 1**

CryLas laser uses one stop bit

**terminator: bytes = b'\n'**

The terminator is LF

**timeout: Union[int, float] = 10**

use 10 seconds timeout as default (a long timeout is needed!)

**class CryLasLaserShutterStatus** (*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)


Bases: `aenum.Enum`

Status of the CryLas laser shutter

**CLOSED = 0**

**OPENED = 1**

## hvl\_ccb.dev.ea\_psi9000

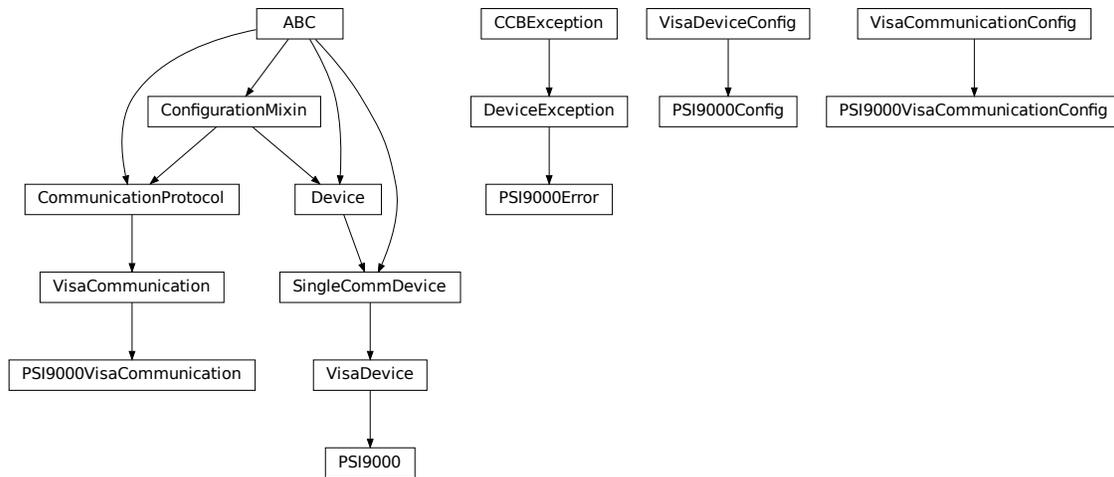

Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to now, all the testing was done with NI-Visa)

```
class PSI9000(com: Union[hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunication,
    hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunicationConfig, dict], dev_config:
    Optional[Union[hvl_ccb.dev.ea_psi9000.PSI9000Config, dict]] = None)
```

Bases: *hvl\_ccb.dev.visa.VisaDevice*

Elektro Automatik PSI 9000 power supply.

**MS\_NOMINAL\_CURRENT** = 2040

**MS\_NOMINAL\_VOLTAGE** = 80

**SHUTDOWN\_CURRENT\_LIMIT** = 0.1

**SHUTDOWN\_VOLTAGE\_LIMIT** = 0.1

**check\_master\_slave\_config()** → *None*

Checks if the master / slave configuration and initializes if successful

**Raises** *PSI9000Error* – if master-slave configuration failed

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns** the VisaCommunication class

**get\_output()** → bool

Reads the current state of the DC output of the source. Returns True, if it is enabled, false otherwise.

**Returns** the state of the DC output

**get\_system\_lock()** → bool

Get the current lock state of the system. The lock state is true, if the remote control is active and false, if not.

**Returns** the current lock state of the device

**get\_ui\_lower\_limits()** → Tuple[float, float]

Get the lower voltage and current limits. A lower power limit does not exist.

**Returns** Umin in V, Imin in A

**get\_ui\_upper\_limits()** → Tuple[float, float, float]

Get the upper voltage, current and power limits.

**Returns** Umax in V, Imax in A, Pmax in W

**get\_voltage\_current\_setpoint()** → Tuple[float, float]

Get the voltage and current setpoint of the current source.

**Returns** Uset in V, Iset in A

**measure\_voltage\_current()** → Tuple[float, float]

Measure the DC output voltage and current

**Returns** Umeas in V, Imeas in A

**set\_lower\_limits**(*voltage\_limit: Optional[float] = None, current\_limit: Optional[float] = None*) → *None*

Set the lower limits for voltage and current. After writing the values a check is performed if the values are set correctly.

**Parameters**

- **voltage\_limit** – is the lower voltage limit in V
- **current\_limit** – is the lower current limit in A

**Raises** *PSI9000Error* – if the limits are out of range

**set\_output**(*target\_onstate: bool*) → *None*

Enables / disables the DC output.

**Parameters** **target\_onstate** – enable or disable the output power

**Raises** *PSI9000Error* – if operation was not successful

**set\_system\_lock**(*lock: bool*) → *None*

Lock / unlock the device, after locking the control is limited to this class unlocking only possible when voltage and current are below the defined limits

**Parameters** **lock** – True: locking, False: unlocking

**set\_upper\_limits**(*voltage\_limit: Optional[float] = None, current\_limit: Optional[float] = None, power\_limit: Optional[float] = None*) → *None*

Set the upper limits for voltage, current and power. After writing the values a check is performed if the values are set. If a parameter is left blank, the maximum configurable limit is set.

**Parameters**

- **voltage\_limit** – is the voltage limit in V
- **current\_limit** – is the current limit in A
- **power\_limit** – is the power limit in W

**Raises** *PSI9000Error* – if limits are out of range

**set\_voltage\_current**(*volt: float, current: float*) → *None*

Set voltage and current setpoints.

After setting voltage and current, a check is performed if writing was successful.

#### Parameters

- **volt** – is the setpoint voltage: 0..81.6 V (1.02 \* 0-80 V) (absolute max, can be smaller if limits are set)
- **current** – is the setpoint current: 0..2080.8 A (1.02 \* 0 - 2040 A) (absolute max, can be smaller if limits are set)

Raises **PSI9000Error** – if the desired setpoint is out of limits

**start**() → *None*

Start this device.

**stop**() → *None*

Stop this device. Turns off output and lock, if enabled.

```
class PSI9000Config(spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2,  
power_limit: Union[int, float] = 43500, voltage_lower_limit: Union[int, float] = 0.0,  
voltage_upper_limit: Union[int, float] = 10.0, current_lower_limit: Union[int, float] = 0.0,  
current_upper_limit: Union[int, float] = 2040.0, wait_sec_system_lock: Union[int, float]  
= 0.5, wait_sec_settings_effect: Union[int, float] = 1, wait_sec_initialisation: Union[int,  
float] = 2)
```

Bases: [hvl\\_ccb.dev.visa.VisaDeviceConfig](#)

Elektro Automatik PSI 9000 power supply device class. The device is communicating over a VISA TCP socket.

Using this power supply, DC voltage and current can be supplied to a load with up to 2040 A and 80 V (using all four available units in parallel). The maximum power is limited by the grid, being at 43.5 kW available through the CEE63 power socket.

**clean\_values**() → *None*

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**current\_lower\_limit: Union[int, float] = 0.0**

Lower current limit in A, depending on the experimental setup.

**current\_upper\_limit: Union[int, float] = 2040.0**

Upper current limit in A, depending on the experimental setup.

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**power\_limit: Union[int, float] = 43500**

Power limit in W depending on the experimental setup. With 3x63A, this is 43.5kW. Do not change this value, if you do not know what you are doing. There is no lower power limit.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**voltage\_lower\_limit: Union[int, float] = 0.0**

Lower voltage limit in V, depending on the experimental setup.

**voltage\_upper\_limit: Union[int, float] = 10.0**

Upper voltage limit in V, depending on the experimental setup.

**wait\_sec\_initialisation: Union[int, float] = 2**

**wait\_sec\_settings\_effect: Union[int, float] = 1**

**wait\_sec\_system\_lock: Union[int, float] = 0.5**

**exception PSI9000Error**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

Base error class regarding problems with the PSI 9000 supply.

**class PSI9000VisaCommunication(configuration)**

Bases: [hvl\\_ccb.comm.visa.VisaCommunication](#)

Communication protocol used with the PSI 9000 power supply.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class PSI9000VisaCommunicationConfig**(*host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], interface\_type: Union[str, hvl\_ccb.comm.visa.VisaCommunicationConfig.InterfaceType] = InterfaceType.TCPIP\_SOCKET, board: int = 0, port: int = 5025, timeout: int = 5000, chunk\_size: int = 204800, open\_timeout: int = 1000, write\_termination: str = '\n', read\_termination: str = '\n', visa\_backend: str = ''*)

Bases: [hvl\\_ccb.comm.visa.VisaCommunicationConfig](#)

Visa communication protocol config dataclass with specification for the PSI 9000 power supply.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign



**property current:** *hvl\_ccb.dev.fug.FuGProbusVSetRegisters*

Returns the registers for the current output

**Returns**

**property current\_monitor:** *hvl\_ccb.dev.fug.FuGProbusVMonitorRegisters*

Returns the registers for the current monitor.

A typically usage will be “self.current\_monitor.value” to measure the output current

**Returns**

**property di:** *hvl\_ccb.dev.fug.FuGProbusVDIRegisters*

Returns the registers for the digital inputs

**Returns** FuGProbusVDIRegisters

**identify\_device()** → *None*

Identify the device nominal voltage and current based on its model number.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**property max\_current:** `Union[int, float]`

Returns the maximal current which could provided within the test setup

**Returns**

**property max\_current\_hardware:** `Union[int, float]`

Returns the maximal current which could provided with the power supply

**Returns**

**property max\_voltage:** `Union[int, float]`

Returns the maximal voltage which could provided within the test setup

**Returns**

**property max\_voltage\_hardware:** `Union[int, float]`

Returns the maximal voltage which could provided with the power supply

**Returns**

**property on:** *hvl\_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the output switch to turn the output on or off

**Returns** FuGProbusVDORegisters

**property outX0:** *hvl\_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X0

**Returns** FuGProbusVDORegisters

**property outX1:** *hvl\_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X1

**Returns** FuGProbusVDORegisters

**property outX2:** *hvl\_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X2

**Returns** FuGProbusVDORegisters

**property outXCMD:** *hvl\_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital outputX-CMD

**Returns** FuGProbusVDORegisters

**start**(*max\_voltage=0, max\_current=0*) → *None*

Opens the communication protocol and configures the device.

**Parameters**

- **max\_voltage** – Configure here the maximal permissible voltage which is allowed in the given experimental setup
- **max\_current** – Configure here the maximal permissible current which is allowed in the given experimental setup

**property voltage:** *hvl\_ccb.dev.fug.FuGProbusVSetRegisters*

Returns the registers for the voltage output

**Returns**

**property voltage\_monitor:** *hvl\_ccb.dev.fug.FuGProbusVMonitorRegisters*

Returns the registers for the voltage monitor.

A typically usage will be “self.voltage\_monitor.value” to measure the output voltage

**Returns**

**class FuGConfig**(*wait\_sec\_stop\_commands: Union[int, float] = 0.5*)

Bases: object

Device configuration dataclass for FuG power supplies.

**clean\_values**()

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**wait\_sec\_stop\_commands: Union[int, float] = 0.5**

Time to wait after subsequent commands during stop (in seconds)

```
class FuGDigitalVal(value)
```

```
    Bases: enum.IntEnum
```

```
    An enumeration.
```

```
    NO = 0
```

```
    OFF = 0
```

```
    ON = 1
```

```
    YES = 1
```

```
exception FuGError(*args, **kwargs)
```

```
    Bases: hvl_ccb.dev.base.DeviceException
```

```
    Error with the FuG voltage source.
```

```
    errorcode: str
```

```
        Errorcode from the Probus, see documentation of Probus V chapter 5. Errors with three-digit errorcodes are thrown by this python module.
```

```
class FuErrorcodes(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

```
    Bases: hvl_ccb.utils.enum.NameEnum
```

```
    The power supply can return an errorcode. These errorcodes are handled by this class. The original errorcodes from the source are with one or two digits, see documentation of Probus V chapter 5. All three-digit errorcodes are from this python module.
```

```
E0 = ('no error', 'standard response on each command')
```

```
E1 = ('no data available', 'Customer tried to read from GPIB but there were no data prepared. (IBIG50 sent command ~T2 to ADDA)')
```

```
E10 = ('unknown SCPI command', 'This SCPI command is not implemented')
```

```
E100 = ('Command is not implemented', 'You tried to execute a command, which is not implemented or does not exist')
```

```
E106 = ('The rampstate is a read-only register', 'You tried to write data to the register, which can only give you the status of the ramping.')
```

```
E11 = ('not allowed Trigger-on-Talk', 'Not allowed attempt to Trigger-on-Talk (~T1) while ADDA was in addressable mode.')
```

```
E115 = ('The given index to select a digital value is out of range', 'Only integer values between 0 and 1 are allowed.')
```

```
E12 = ('invalid argument in ~Tn command', 'Only ~T1 and ~T2 is implemented.')
```

```
E125 = ('The given index to select a ramp mode is out of range', 'Only integer values between 0 and 4 are allowed.')
```

```
E13 = ('invalid N-value', 'Register > K8 contained an invalid value. Error code is output on an attempt to query data with ? or ~T1')
```

```
E135 = ('The given index to select the readback channel is out of range', 'Only integer values between 0 and 6 are allowed.')
```

```
E14 = ('register is write only', 'Some registers can only be writte to (i.e.> H0)')
```

```
E145 = ('The given value for the AD-conversion is unknown', 'Valid values for the ad-conversion are integer values from "0" to "7".')
```

```
E15 = ('string too long', 'i.e.serial number string too long during calibration')
```

E155 = ('The given value to select a polarity is out range.', 'The value should be 0 or 1.')  
E16 = ('wrong checksum', 'checksum over command string was not correct, refer also to 4.4 of the Probus V documentation')  
E165 = ('The given index to select the terminator string is out of range', '')  
E2 = ('unknown register type', "No valid register type after '>'")  
E206 = ('This status register is read-only', 'You tried to write data to this register, which can only give you the actual status of the corresponding digital output.')  
E306 = ('The monitor register is read-only', 'You tried to write data to a monitor, which can only give you measured data.')  
E4 = ('invalid argument', 'The argument of the command was rejected .i.e. malformed number')  
E5 = ('argument out of range', 'i.e. setvalue higher than type value')  
E504 = ('Empty string as response', 'The connection is broken.')  
E505 = ('The returned register is not the requested.', 'Maybe the connection is overburden.')  
E6 = ('register is read only', 'Some registers can only be read but not written to. (i.e. monitor registers)')  
E666 = ('You cannot overwrite the most recent error in the interface of the power supply. But, well: You created an error anyway...', '')  
E7 = ('Receive Overflow', 'Command string was longer than 50 characters.')  
E8 = ('EEPROM is write protected', 'Write attempt to calibration data while the write protection switch was set to write protected.')  
E9 = ('address error', 'A non addressed command was sent to ADDA while it was in addressable mode (and vice versa).')  
raise\_()

**class** FuGMonitorModes(*value*)

Bases: enum.IntEnum

An enumeration.

**T1MS** = 1

15 bit + sign, 1 ms integration time

**T200MS** = 6

typ. 19 bit + sign, 200 ms integration time

**T20MS** = 3

17 bit + sign, 20 ms integration time

**T256US** = 0

14 bit + sign, 256 us integration time

**T40MS** = 4

17 bit + sign, 40 ms integration time

**T4MS** = 2

15 bit + sign, 4 ms integration time

**T800MS = 7**  
typ. 20 bit + sign, 800 ms integration time

**T80MS = 5**  
typ. 18 bit + sign, 80 ms integration time

**class FuGPolarities**(*value*)

Bases: `enum.IntEnum`

An enumeration.

**NEGATIVE = 1**

**POSITIVE = 0**

**class FuGProbusIV**(*com*, *dev\_config=None*)

Bases: `hvl_ccb.dev.base.SingleCommDevice`, `abc.ABC`

FuG Probus IV device class

Sends basic SCPI commands and reads the answer. Only the special commands and PROBUS IV instruction set is implemented.

**command**(*command*: `hvl_ccb.dev.fug.FuGProbusIVCommands`, *value=None*) → str

#### Parameters

- **command** – one of the commands given within `FuGProbusIVCommands`
- **value** – an optional value, depending on the command

**Returns** a String if a query was performed

**static config\_cls**()

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls**()

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**output\_off**() → *None*

Switch DC voltage output off.

**reset**() → *None*

Reset of the interface: All setvalues are set to zero

**abstract start**()

Open the associated communication protocol.

**stop**() → *None*

Close the associated communication protocol.

**class FuGProbusIVCommands**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.NameEnum`

An enumeration.

**ADMODE = ('S', (<enum 'FuGMonitorModes'>, <class 'int'>))**

**CURRENT = ('I', (<class 'int'>, <class 'float'>))**

**EXECUTE = ('X', None)**

**EXECUTEONX** = ('G', (<enum 'FuGDigitalVal'>, <class 'int'>))

Wait for “X” to execute pending commands

**ID** = (\*IDN?, None)

**OUTPUT** = ('F', (<enum 'FuGDigitalVal'>, <class 'int'>))

**POLARITY** = ('P', (<enum 'FuGPolarities'>, <class 'int'>))

**QUERY** = ('?', None)

**READBACKCHANNEL** = ('N', (<enum 'FuGReadbackChannels'>, <class 'int'>))

**RESET** = ('=', None)

**TERMINATOR** = ('Y', (<enum 'FuGTerminators'>, <class 'int'>))

**VOLTAGE** = ('U', (<class 'int'>, <class 'float'>))

**XOUTPUTS** = ('R', <class 'int'>)

TODO: the possible values are limited to 0..13

**class FuGProbusV**(*com, dev\_config=None*)

Bases: [hvl\\_ccb.dev.fug.FuGProbusIV](#)

FuG Probus V class which uses register based commands to control the power supplies

**get\_register**(*register: str*) → str

get the value from a register

**Parameters** **register** – the register from which the value is requested

**Returns** the value of the register as a String

**set\_register**(*register: str, value: Union[int, float, str]*) → None

generic method to set value to register

**Parameters**

- **register** – the name of the register to set the value
- **value** – which should be written to the register

**class FuGProbusVConfigRegisters**(*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Configuration and Status values, acc. 4.2.5

**property execute\_on\_x:** [hvl\\_ccb.dev.fug.FuGDigitalVal](#)

status of Execute-on-X

**Returns** FuGDigitalVal of the status

**property most\_recent\_error:** [hvl\\_ccb.dev.fug.FuErrorcodes](#)

Reads the Error-Code of the most recent command

**Return** FuGError

**Raises** [FuGError](#) – if code is not “E0”

**property readback\_data:** [hvl\\_ccb.dev.fug.FuGReadbackChannels](#)

Preselection of readout data for Trigger-on-Talk

**Returns** index for the readback channel

**property srq\_mask:** int

SRQ-Mask, Service-Request Enable status bits for SRQ 0: no SRQ Bit 2: SRQ on change of status to CC  
Bit 1: SRQ on change to CV

**Returns** representative integer value

**property srq\_status:** **str**

SRQ-Statusbyte output as a decimal number: Bit 2: PS is in CC mode Bit 1: PS is in CV mode

**Returns** representative string

**property status:** **str**

Statusbyte as a string of 0/1. Combined status (compatibel to Probus IV), MSB first: Bit 7: I-REG Bit 6: V-REG Bit 5: ON-Status Bit 4: 3-Reg Bit 3: X-Stat (polarity) Bit 2: Cal-Mode Bit 1: unused Bit 0: SEL-D

**Returns** string of 0/1

**property terminator:** *hvl\_ccb.dev.fug.FuGTerminators*

Terminator character for answer strings from ADDA

**Returns** FuGTerminators

**class FuGProbusVDIRegisters**(*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Digital Inputs acc. 4.2.4

**property analog\_control:** *hvl\_ccb.dev.fug.FuGDigitalVal*

**Returns** shows 1 if power supply is controlled by the analog interface

**property calibration\_mode:** *hvl\_ccb.dev.fug.FuGDigitalVal*

**Returns** shows 1 if power supply is in calibration mode

**property cc\_mode:** *hvl\_ccb.dev.fug.FuGDigitalVal*

**Returns** shows 1 if power supply is in CC mode

**property cv\_mode:** *hvl\_ccb.dev.fug.FuGDigitalVal*

**Returns** shows 1 if power supply is in CV mode

**property digital\_control:** *hvl\_ccb.dev.fug.FuGDigitalVal*

**Returns** shows 1 if power supply is digitally controlled

**property on:** *hvl\_ccb.dev.fug.FuGDigitalVal*

**Returns** shows 1 if power supply ON

**property reg\_3:** *hvl\_ccb.dev.fug.FuGDigitalVal*

For special applications.

**Returns** input from bit 3-REG

**property x\_stat:** *hvl\_ccb.dev.fug.FuGPolarities*

**Returns** polarity of HVPS with polarity reversal

**class FuGProbusVDORegisters**(*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Digital outputs acc. 4.2.2

**property out:** Union[int, *hvl\_ccb.dev.fug.FuGDigitalVal*]

Status of the output according to the last setting. This can differ from the actual state if output should only pulse.

**Returns** FuGDigitalVal

**property status:** *hvl\_ccb.dev.fug.FuGDigitalVal*

Returns the actual value of output. This can differ from the set value if pulse function is used.

**Returns** FuGDigitalVal

**class FuGProbusVMonitorRegisters**(*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Analog monitors acc. 4.2.3

**property adc\_mode:** *hvl\_ccb.dev.fug.FuGMonitorModes*

The programmed resolution and integration time of the AD converter

**Returns** FuGMonitorModes

**property value:** float

Value from the monitor.

**Returns** a float value in V or A

**property value\_raw:** float

uncalibrated raw value from AD converter

**Returns** float value from ADC

**class FuGProbusVRegisterGroups**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl\_ccb.utils.enum.NameEnum*

An enumeration.

CONFIG = 'K'

INPUT = 'D'

MONITOR\_I = 'M1'

MONITOR\_V = 'M0'

OUTPUTONCMD = 'BON'

OUTPUTX0 = 'B0'

OUTPUTX1 = 'B1'

OUTPUTX2 = 'B2'

OUTPUTXCMD = 'BX'

SETCURRENT = 'S1'

SETVOLTAGE = 'S0'

**class FuGProbusVSetRegisters**(*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Setvalue control acc. 4.2.1 for the voltage and the current output

**property actualsetvalue:** float

The actual valid set value, which depends on the ramp function.

**Returns** actual valid set value

**property high\_resolution:** *hvl\_ccb.dev.fug.FuGDigitalVal*

Status of the high resolution mode of the output.

**Return 0** normal operation

**Return 1** High Res. Mode

**property rampmode:** *hvl\_ccb.dev.fug.FuGRampModes*

The set ramp mode to control the setvalue.

**Returns** the mode of the ramp as instance of FuGRampModes

**property ramprate:** **float**

The set ramp rate in V/s.

**Returns** ramp rate in V/s

**property rampstate:** *hvl\_ccb.dev.fug.FuGDigitalVal*

Status of ramp function.

**Return 0** if final setvalue is reached

**Return 1** if still ramping up

**property setvalue:** **float**

For the voltage or current output this setvalue was programmed.

**Returns** the programmed setvalue

**class FuGRampModes**(*value*)

Bases: `enum.IntEnum`

An enumeration.

**FOLLOWRAMP = 1**

Follow the ramp up- and downwards

**IMMEDIATELY = 0**

Standard mode: no ramp

**ONLYUPWARDSOFFTOZERO = 4**

Follow the ramp up- and downwards, if output is OFF set value is zero

**RAMPUPWARDS = 2**

Follow the ramp only upwards, downwards immediately

**SPECIALRAMPUPWARDS = 3**

Follow a special ramp function only upwards

**class FuGReadbackChannels**(*value*)

Bases: `enum.IntEnum`

An enumeration.

**CURRENT = 1**

**FIRMWARE = 5**

**RATEDCURRENT = 4**

**RATEDVOLTAGE = 3**

**SN = 6**

**STATUSBYTE = 2**

**VOLTAGE = 0**

**class** `FuGSerialCommunication`(*configuration*)

Bases: `hvl_ccb.comm.serial.SerialCommunication`

Specific communication protocol implementation for FuG power supplies. Already predefines device-specific protocol parameters in config.

**static** `config_cls()`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query**(*command: str*) → str

Send a command to the interface and handle the status message. Eventually raises an exception.

**Parameters** `command` – Command to send

**Raises** `FuGError` – if the connection is broken or the error from the power source itself

**Returns** Answer from the interface or empty string

**class** `FuGSerialCommunicationConfig`(*terminator: bytes = b'\n', encoding: str = 'utf-8', encoding\_error\_handling: str = 'strict', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)*

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate: int = 9600**

Baudrate for FuG power supplies is 9600 baud

**bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8**

One byte is eight bits long

**default\_n\_attempts\_read\_text\_nonempty: int = 10**

default number of attempts to read a non-empty text

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod** `keys()` → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod** `optional_defaults()` → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** Union[str, *hvl\_ccb.comm.serial.SerialCommunicationParity*] = 'N'

FuG does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** Union[int, *hvl\_ccb.comm.serial.SerialCommunicationStopbits*] = 1

FuG uses one stop bit

**terminator:** bytes = b'\n'

The terminator is LF

**timeout:** Union[int, float] = 3

use 3 seconds timeout as default

**wait\_sec\_read\_text\_nonempty:** Union[int, float] = 0.5

default time to wait between attempts of reading a non-empty text

**class FuGTerminators**(*value*)

Bases: enum.IntEnum

An enumeration.


**CR** = 3

**CRLF** = 0

**LF** = 2

**LFCR** = 1

## hvl\_ccb.dev.heinzinger

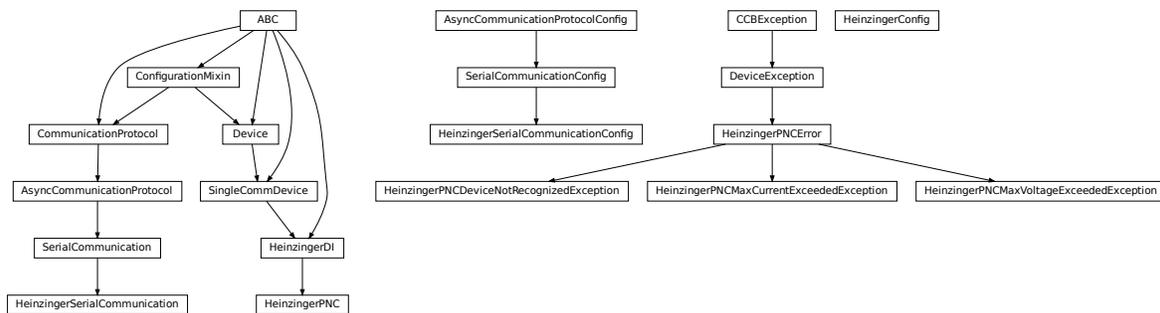

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Manufacturer homepage: <https://www.heinzinger.com/products/accessories-and-more/digital-interfaces/>

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class HeinzingerPNC is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCcap. Manufacturer homepage: <https://www.heinzinger.com/products/high-voltage/universal-high-voltage-power-supplies/>

```
class HeinzingerConfig(default_number_of_recordings: Union[int,  
                       hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum] = 1,  
                       number_of_decimals: int = 6, wait_sec_stop_commands: Union[int, float] = 0.5)
```

Bases: object

Device configuration dataclass for Heinzinger power supplies.

```
class RecordingsEnum(value)
```

Bases: enum.IntEnum

An enumeration.

**EIGHT = 8**

**FOUR = 4**

**ONE = 1**

**SIXTEEN = 16**

**TWO = 2**

```
clean_values()
```

```
default_number_of_recordings: Union[int,  
hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum] = 1
```

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
is_configdataclass = True
```

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

```
number_of_decimals: int = 6
```

```
classmethod optional_defaults() → Dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

```
wait_sec_stop_commands: Union[int, float] = 0.5
```

Time to wait after subsequent commands during stop (in seconds)

```
class HeinzingerDI(com, dev_config=None)
```

Bases: *hvl\_ccb.dev.base.SingleCommDevice*, abc.ABC

Heinzinger Digital Interface I/II device class

Sends basic SCPI commands and reads the answer. Only the standard instruction set from the manual is implemented.

**class OutputStatus**(*value*)

Bases: `enum.IntEnum`

Status of the voltage output

**OFF** = 0

**ON** = 1

**UNKNOWN** = -1

**static config\_cls**()

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls**()

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**get\_current**() → float

Queries the set current of the Heinzinger PNC (not the measured current!).

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**get\_interface\_version**() → str

Queries the version number of the digital interface.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**get\_number\_of\_recordings**() → int

Queries the number of recordings the device is using for average value calculation.

**Returns** int number of recordings

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**get\_serial\_number**() → str

Ask the device for its serial number and returns the answer as a string.

**Returns** string containing the device serial number

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**get\_voltage**() → float

Queries the set voltage of the Heinzinger PNC (not the measured voltage!).

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**measure\_current**() → float

Ask the Device to measure its output current and return the measurement result.

**Returns** measured current as float

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**measure\_voltage**() → float

Ask the Device to measure its output voltage and return the measurement result.

**Returns** measured voltage as float

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**output\_off()** → *None*

Switch DC voltage output off and updates the output status.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**output\_on()** → *None*

Switch DC voltage output on and updates the output status.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**property output\_status:** *hvl\_ccb.dev.heinzinger.HeinzingerDI.OutputStatus*

**reset\_interface()** → *None*

Reset of the digital interface; only Digital Interface I: Power supply is switched to the Local-Mode (Manual operation)

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_current(value: Union[int, float])** → *None*

Sets the output current of the Heinzinger PNC to the given value.

**Parameters value** – current expressed in *self.unit\_current*

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_number\_of\_recordings(value: Union[int, hvl\_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum])** → *None*

Sets the number of recordings the device is using for average value calculation. The possible values are 1, 2, 4, 8 and 16.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_voltage(value: Union[int, float])** → *None*

Sets the output voltage of the Heinzinger PNC to the given value.

**Parameters value** – voltage expressed in *self.unit\_voltage*

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**abstract start()**

Opens the communication protocol.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened.

**stop()** → *None*

Stop the device. Closes also the communication protocol.

**class HeinzingerPNC(com, dev\_config=None)**

Bases: *hvl\_ccb.dev.heinzinger.HeinzingerDI*

Heinzinger PNC power supply device class.

The power supply is controlled over a Heinzinger Digital Interface I/II

**class UnitCurrent(value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None)**

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

An enumeration.

**A** = 3

**UNKNOWN** = 1

**mA** = 2

```

class UnitVoltage(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl_ccb.utils.enum.AutoNumberNameEnum

    An enumeration.

    UNKNOWN = 1

    V = 2

    kV = 3

identify_device() → None
    Identify the device nominal voltage and current based on its serial number.

    Raises SerialCommunicationIOError – when communication port is not opened

property max_current: Union[int, float]
property max_current_hardware: Union[int, float]
property max_voltage: Union[int, float]
property max_voltage_hardware: Union[int, float]
set_current(value: Union[int, float]) → None
    Sets the output current of the Heinzinger PNC to the given value.

    Parameters value – current expressed in self.unit_current

    Raises SerialCommunicationIOError – when communication port is not opened

set_voltage(value: Union[int, float]) → None
    Sets the output voltage of the Heinzinger PNC to the given value.

    Parameters value – voltage expressed in self.unit_voltage

    Raises SerialCommunicationIOError – when communication port is not opened

start() → None
    Opens the communication protocol and configures the device.

property unit_current: hvl_ccb.dev.heinzinger.HeinzingerPNC.UnitCurrent
property unit_voltage: hvl_ccb.dev.heinzinger.HeinzingerPNC.UnitVoltage

exception HeinzingerPNCDeviceNotRecognizedException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

    Error indicating that the serial number of the device is not recognized.

exception HeinzingerPNCError
    Bases: hvl_ccb.dev.base.DeviceException

    General error with the Heinzinger PNC voltage source.

exception HeinzingerPNCMaxCurrentExceededException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

    Error indicating that program attempted to set the current to a value exceeding ‘max_current’.

exception HeinzingerPNCMaxVoltageExceededException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

    Error indicating that program attempted to set the voltage to a value exceeding ‘max_voltage’.

class HeinzingerSerialCommunication(configuration)
    Bases: hvl_ccb.comm.serial.SerialCommunication

```

Specific communication protocol implementation for Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class HeinzingerSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 40, port:
                                         Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                         Union[str, hvl_ccb.comm.serial.SerialCommunicationParity]
                                         = <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                         Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                         <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                         Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                         <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                         Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate: int = 9600**

Baudrate for Heinzinger power supplies is 9600 baud

**bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8**

One byte is eight bits long

**default\_n\_attempts\_read\_text\_nonempty: int = 40**

increased to 40 default number of attempts to read a non-empty text

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'**

Heinzinger does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


**stopbits:** `Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1`  
Heinzinger uses one stop bit

**terminator:** `bytes = b'\n'`  
The terminator is LF

**timeout:** `Union[int, float] = 3`  
use 3 seconds timeout as default

**wait\_sec\_read\_text\_nonempty:** `Union[int, float] = 0.5`  
default time to wait between attempts of reading a non-empty text

### `hvl_ccb.dev.labjack`

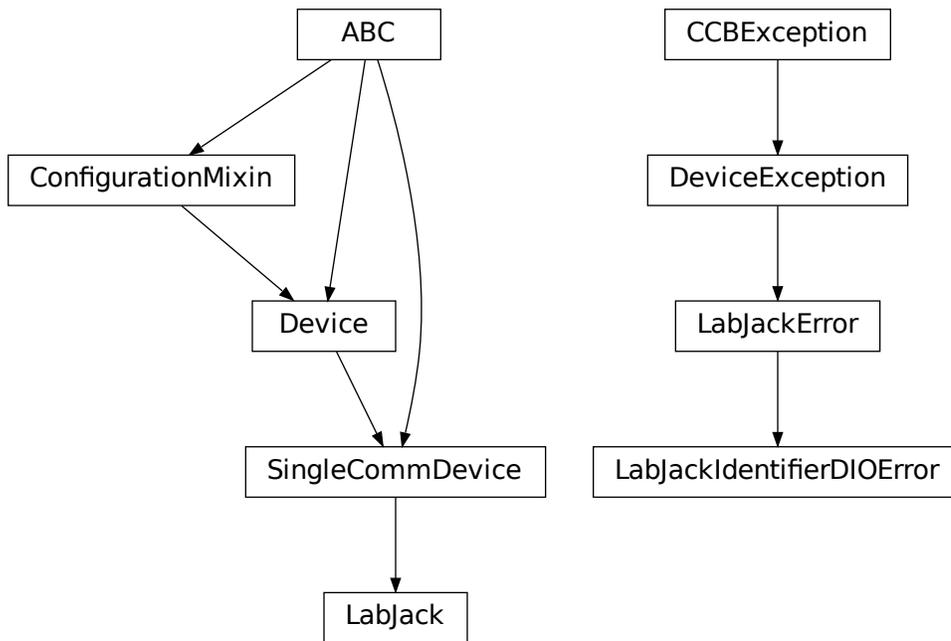

LabJack T-series devices wrapper around the LabJack's LJM Library; see <https://labjack.com/ljm> . The wrapper was originally developed and tested for a LabJack T7-PRO device. A

## Extra installation

To use this LabJack T-series devices wrapper:

1. install the `hvl_ccb` package with a `labjack` extra feature:

```
$ pip install "hvl_ccb[labjack]"
```

this will install the Python bindings for the library.

2. install the library - follow instruction in <https://labjack.com/support/software/installers/ljm> .

**class LabJack**(*com, dev\_config=None*)

Bases: `hvl_ccb.dev.base.SingleCommDevice`

LabJack Device.

This class is tested with a LabJack T7-Pro and should also work with T4 and T7 devices communicating through the LJM Library. Other or older hardware versions and variants of LabJack devices are not supported.

**class AInRange**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `hvl_ccb.utils.enum.StrEnumBase`

An enumeration.

**ONE = 1.0**

**ONE\_HUNDREDTH = 0.01**

**ONE\_TENTH = 0.1**

**TEN = 10.0**

**property value: float**

**class BitLimit**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `aenum.IntEnum`

Maximum integer values for clock settings

**THIRTY\_TWO\_BIT = 4294967295**

**class CalMicroAmpere**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `aenum.Enum`

Pre-defined microampere (uA) values for calibration current source query.

**TEN = '10uA'**

**TWO\_HUNDRED = '200uA'**

**class CjcType**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `hvl_ccb.utils.enum.NameEnum`

CJC slope and offset

**internal = (1, 0)**

**lm34 = (55.56, 255.37)**

**class ClockFrequency**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `aenum.IntEnum`

Available clock frequencies, in Hz

```

FIVE_MHZ = 5000000
FORTY_MHZ = 40000000
MAXIMUM = 80000000
MINIMUM = 312500
TEN_MHZ = 10000000
TWELVE_HUNDRED_FIFTY_KHZ = 1250000
TWENTY_FIVE_HUNDRED_KHZ = 2500000
TWENTY_MHZ = 20000000

```

**DIOChannel**

alias of `hvl_ccb._dev.labjack.TSeriesDIOChannel`

```
class DIOStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `aenum.IntEnum`

State of a digital I/O channel.

```
HIGH = 1
```

```
LOW = 0
```

```
class DeviceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `hvl_ccb.utils.enum.AutoNumberNameEnum`

LabJack device types.

Can be also looked up by ambiguous Product ID (*p\_id*) or by instance name: ``python LabJackDeviceType(4) is LabJackDeviceType('T4')``

```
ANY = 1
```

```
T4 = 2
```

```
T7 = 3
```

```
T7_PRO = 4
```

```
classmethod get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
List[hvl_ccb._dev.labjack.DeviceType]]
```

Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

**Parameters** *p\_id* – Product ID of a LabJack device

**Returns** Instance or list of instances of `LabJackDeviceType`

**Raises** `ValueError` – when Product ID is unknown

```
class TemperatureUnit(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
```

Bases: `hvl_ccb.utils.enum.NameEnum`

Temperature unit (to be returned)

```
C = 1
```

```
F = 2
```

```
K = 0
```

```
class ThermocoupleType(value=<no_arg>, names=None, module=None, type=None, start=1,  
                        boundary=None)
```

Bases: `hvl_ccb.utils.enum.NameEnum`

Thermocouple type; NONE means disable thermocouple mode.

**C = 30**

**E = 20**

**J = 21**

**K = 22**

**NONE = 0**

**PT100 = 40**

**PT1000 = 42**

**PT500 = 41**

**R = 23**

**S = 25**

**T = 24**

```
config_high_pulse(address: Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel], t_start: Union[int,  
float], t_width: Union[int, float], n_pulses: int = 1) → None
```

Configures one FIO channel to send a timed HIGH pulse. Configure multiple channels to send pulses with relative timing accuracy. Times have a maximum resolution of 1e-7 seconds @ 10 MHz. :param address: FIO channel: [T7] FIO0;2;3;4;5. [T4] FIO6;7. :raises LabJackError if address is not supported. :param t\_start: pulse start time in seconds. :raises ValueError: if t\_start is negative or would exceed the clock period. :param t\_width: duration of high pulse, in seconds. :raises ValueError: if t\_width is negative or would exceed the clock period. :param n\_pulses: number of pulses to be sent; single pulse default. :raises TypeError if n\_pulses is not of type int. :raises Value Error if n\_pulses is negative or exceeds the 32bit limit.

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

```
disable_pulses(*addresses: Optional[Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel]]) → None
```

Disable previously configured pulse channels. :param addresses: tuple of FIO addresses. All channels disabled if no argument is passed.

```
enable_clock(clock_enabled: bool) → None
```

Enable/disable LabJack clock to configure or send pulses. :param clock\_enabled: True -> enable, False -> disable. :raises TypeError: if clock\_enabled is not of type bool

```
get_ain(*channels: int) → Union[float, Sequence[float]]
```

Read currently measured value (voltage, resistance, ...) from one or more of analog inputs.

**Parameters channels** – AIN number or numbers (0..254)

**Returns** the read value (voltage, resistance, ...) as *float* or *tuple* of them in case multiple channels given

```
get_cal_current_source(name: Union[str, CalMicroAmpere]) → float
```

This function will return the calibration of the chosen current source, this is not a measurement!

The value was stored during fabrication.

**Parameters** **name** – ‘200uA’ or ‘10uA’ current source

**Returns** calibration of the chosen current source in ampere

**get\_clock()** → Dict[str, Union[int, float]]

Return clock settings read from LabJack.

**get\_digital\_input**(*address: Union[str, hvl\_ccb.\_dev.labjack.TSeriesDIOChannel]*) → *hvl\_ccb.dev.labjack.LabJack.DIOStatus*

Get the value of a digital input.

allowed names for T7 (Pro): FIO0 - FIO7, EIO0 - EIO 7, CIO0- CIO3, MIO0 - MIO2 :param address: name of the output -> ‘FIO0’ :return: HIGH when *address* DIO is high, and LOW when *address* DIO is low

**get\_product\_id()** → int

This function returns the product ID reported by the connected device.

Attention: returns 7 for both T7 and T7-Pro devices!

**Returns** integer product ID of the device

**get\_product\_name**(*force\_query\_id=False*) → str

This function will return the product name based on product ID reported by the device.

Attention: returns “T7” for both T7 and T7-Pro devices!

**Parameters** **force\_query\_id** – boolean flag to force *get\_product\_id* query to device instead of using cached device type from previous queries.

**Returns** device name string, compatible with *LabJack.DeviceType*

**get\_product\_type**(*force\_query\_id: bool = False*) → *hvl\_ccb.\_dev.labjack.DeviceType*

This function will return the device type based on reported device type and in case of unambiguity based on configuration of device’s communication protocol (e.g. for “T7” and “T7\_PRO” devices), or, if not available first matching.

**Parameters** **force\_query\_id** – boolean flag to force *get\_product\_id* query to device instead of using cached device type from previous queries.

**Returns** *DeviceType* instance

**Raises** *LabJackIdentifierDIOError* – when read Product ID is unknown

**get\_sbush\_rh**(*number: int*) → float

Read the relative humidity value from a serial SBUS sensor.

**Parameters** **number** – port number (0..22)

**Returns** relative humidity in %RH

**get\_sbush\_temp**(*number: int*) → float

Read the temperature value from a serial SBUS sensor.

**Parameters** **number** – port number (0..22)

**Returns** temperature in Kelvin

**get\_serial\_number()** → int

Returns the serial number of the connected LabJack.

**Returns** Serial number.

**read\_resistance**(*channel: int*) → float

Read resistance from specified channel.

**Parameters** `channel` – channel with resistor

**Returns** resistance value with 2 decimal places

**read\_thermocouple**(*pos\_channel: int*) → float

Read the temperature of a connected thermocouple.

**Parameters** `pos_channel` – is the AIN number of the positive pin

**Returns** temperature in specified unit

**send\_pulses**(*\*addresses: Union[str, hvl\_ccb.\_dev.labjack.TSeriesDIOChannel]*) → None

Sends pre-configured pulses for specified addresses. :param addresses: tuple of FIO addresses :raises LabJackError if an address has not been configured.

**set\_ain\_differential**(*pos\_channel: int, differential: bool*) → None

Sets an analog input to differential mode or not. T7-specific: For base differential channels, positive must be even channel from 0-12 and negative must be positive+1. For extended channels 16-127, see Mux80 datasheet.

**Parameters**

- `pos_channel` – is the AIN number (0..12)
- `differential` – True or False

**Raises** `LabJackError` – if parameters are unsupported

**set\_ain\_range**(*channel: int, vrange: Union[Real, AInRange]*) → None

Set the range of an analog input port.

**Parameters**

- `channel` – is the AIN number (0..254)
- `vrange` – is the voltage range to be set

**set\_ain\_resistance**(*channel: int, vrange: Union[Real, AInRange], resolution: int*) → None

Set the specified channel to resistance mode. It utilized the 200uA current source of the LabJack.

**Parameters**

- `channel` – channel that should measure the resistance
- `vrange` – voltage range of the channel
- `resolution` – resolution index of the channel T4: 0-5, T7: 0-8, T7-Pro 0-12

**set\_ain\_resolution**(*channel: int, resolution: int*) → None

Set the resolution index of an analog input port.

**Parameters**

- `channel` – is the AIN number (0..254)
- `resolution` – is the resolution index within 0...`get\_product\_type().ain\_max\_resolution` range; 0 will set the resolution index to default value.

**set\_ain\_thermocouple**(*pos\_channel: int, thermocouple: Union[None, str, ThermocoupleType], cjc\_address: int = 60050, cjc\_type: Union[str, CjcType] = CjcType.internal, vrange: Union[Real, AInRange] = AInRange.ONE\_HUNDREDTH, resolution: int = 10, unit: Union[str, TemperatureUnit] = TemperatureUnit.K*) → None

Set the analog input channel to thermocouple mode.

**Parameters**

- `pos_channel` – is the analog input channel of the positive part of the differential pair

- **thermocouple** – None to disable thermocouple mode, or string specifying the thermocouple type
- **cjc\_address** – modbus register address to read the CJC temperature
- **cjc\_type** – determines cjc slope and offset, ‘internal’ or ‘lm34’
- **vrange** – measurement voltage range
- **resolution** – resolution index (T7-Pro: 0-12)
- **unit** – is the temperature unit to be returned (‘K’, ‘C’ or ‘F’)

Raises **LabJackError** – if parameters are unsupported

**set\_analog\_output**(*channel: int, value: Union[int, float]*) → *None*

Set the voltage of an analog output port

#### Parameters

- **channel** – DAC channel number 1/0
- **value** – The output voltage value 0-5 Volts int/float

**set\_clock**(*clock\_frequency: Union[Number, ClockFrequency] = 10000000, clock\_period: Number = 1*) → *None*

Configure LabJack clock for pulse out feature. :param clock\_frequency: clock frequency in Hz; default 10 MHz for base 10. :raises ValueError: if clock\_frequency is not allowed (see ClockFrequency). :param clock\_period: clock roll time in seconds; default 1s, 0 for max. :raises ValueError: if clock\_period exceeds the 32bit tick limit. Clock period determines pulse spacing when using multi-pulse settings. Ensure period exceeds maximum intended pulse end time.

**set\_digital\_output**(*address: str, state: Union[int, DIOStatus]*) → *None*

Set the value of a digital output.

#### Parameters

- **address** – name of the output -> ‘FIO0’
- **state** – state of the output -> *DIOStatus* instance or corresponding *int* value

**start**() → *None*

Start the Device.

**stop**() → *None*

Stop the Device.

**exception LabJackError**

Bases: *hvl\_ccb.dev.base.DeviceException*


General Exception for the LabJack device.

**exception LabJackIdentifierDIOError**

Bases: *hvl\_ccb.dev.labjack.LabJackError*

Error indicating a wrong DIO identifier

**hvl\_ccb.dev.lauda**

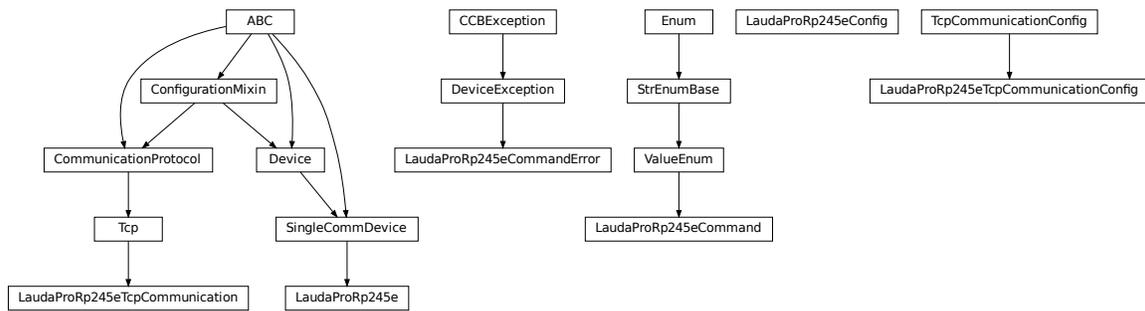

Device class for controlling a Lauda PRO RP245E, circulation chiller over TCP.

**class** `LaudaProRp245e`(*com, dev\_config=None*)

Bases: `hvl_ccb.dev.base.SingleCommDevice`

Lauda RP245E circulation chiller class.

**static config\_cls**() → `Type[hvl_ccb.dev.lauda.LaudaProRp245eConfig]`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**continue\_ramp**() → str

Continue current ramp program.

**Returns** reply of the device to the last call of “query”

**static default\_com\_cls**() → `Type[hvl_ccb.dev.lauda.LaudaProRp245eTcpCommunication]`

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**get\_bath\_temp**() → float

:return : float value of measured lauda bath temp in °C

**get\_device\_type**() → str

:return : Connected Lauda device type (for connection/com test)

**pause**() → str

Stop temperature control and pump.

**Returns** reply of the device to the last call of “query”

**pause\_ramp**() → str

Pause current ramp program.

**Returns** reply of the device to the last call of “query”

**reset\_ramp**() → str

Delete all segments from current ramp program.

**Returns** reply of the device to the last call of “query”

**run**() → str

Start temperature control & pump.

**Returns** reply of the device to the last call of “query”

**set\_control\_mode**(*mod: Union[int, hvl\_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = ExtControlModeEnum.INTERNAL*) → str

Define control mode. 0 = INTERNAL (control bath temp), 1 = EXPT100 (pt100 attached to chiller), 2 = ANALOG, 3 = SERIAL, 4 = USB, 5 = ETH (to be used when passing the ext. temp. via ethernet) (temperature then needs to be passed every second, when not using options 3, 4, or 5)

**Parameters** **mod** – temp control mode (control internal temp or external temp).

**Returns** reply of the device to the last call of “query” (“OK”, if command was recognized”)

**set\_external\_temp**(*external\_temp: float = 20.0*) → str

Pass value of external controlled temperature. Should be done every second, when control of external temperature is active. Has to be done right before control of external temperature is activated.

**Parameters** **external\_temp** – current value of external temperature to be controlled.

**Returns** reply of the device to the last call of “query”

**set\_pump\_level**(*pump\_level: int = 6*) → str

Set pump level Raises ValueError, if pump level is invalid.

**Parameters** **pump\_level** – pump level.

**Returns** reply of the device to the last call of “query”

**set\_ramp\_iterations**(*num: int = 1*) → str

Define number of ramp program cycles.

**Parameters** **num** – number of program cycles to be performed.

**Returns** reply of the device to the last call of “query”

**set\_ramp\_program**(*program: int = 1*) → str

Define ramp program for following ramp commands. Raises ValueError if maximum number of ramp programs (5) is exceeded.

**Parameters** **program** – Number of ramp program to be activated for following commands.

**Returns** reply of the device to the last call of “query”

**set\_ramp\_segment**(*temp: float = 20.0, dur: int = 0, tol: float = 0.0, pump: int = 6*) → str

Define segment of current ramp program - will be attached to current program. Raises ValueError, if pump level is invalid.

**Parameters**

- **temp** – target temperature of current ramp segment
- **dur** – duration in minutes, in which target temperature should be reached
- **tol** – tolerance at which target temperature should be reached (for 0.00, next segment is started after dur has passed).
- **pump** – pump level to be used for this program segment.

**Returns** reply of the device to the last call of “query”

**set\_temp\_set\_point**(*temp\_set\_point: float = 20.0*) → str

Define temperature set point

**Parameters** **temp\_set\_point** – temperature set point.

**Returns** reply of the device to the last call of “query”

**start()** → *None*

Start this device.

**start\_ramp()** → str

Start current ramp program.

**Returns** reply of the device to the last call of “query”

**stop()** → *None*

Stop this device. Disables access and closes the communication protocol.

**stop\_ramp()** → str

Stop current ramp program.

**Returns** reply of the device to the last call of “query”

**validate\_pump\_level**(*level: int*)

Validates pump level. Raises ValueError, if pump level is incorrect. :param level: pump level, integer

**class** **LaudaProRp245eCommand**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl\_ccb.utils.enum.ValueEnum*

Commands for Lauda PRO RP245E Chiller Command strings most often need to be complimented with a parameter (attached as a string) before being sent to the device. Commands implemented as defined in “Lauda Betriebsanleitung fuer PRO Badthermostate und Umwaelzthermostate” pages 42 - 49

**BATH\_TEMP** = 'IN\_PV\_00'

Request internal bath temperature

**COM\_TIME\_OUT** = 'OUT\_SP\_08\_'

Define communication time out

**CONT\_MODE** = 'OUT\_MODE\_01\_'

Set control mode 1=internal, 2=ext. analog, 3=ext. serial, 4=USB, 5=ethernet

**DEVICE\_TYPE** = 'TYPE'

Request device type

**EXTERNAL\_TEMP** = 'OUT\_PV\_05\_'

Pass on external controlled temperature

**LOWER\_TEMP** = 'OUT\_SP\_05\_'

Define lower temp limit

**OPERATION\_MODE** = 'OUT\_SP\_02\_'

Define operation mode

**PUMP\_LEVEL** = 'OUT\_SP\_01\_'

Define pump level 1-8

**RAMP\_CONTINUE** = 'RMP\_CONT'

Continue a paused ramp program

**RAMP\_DELETE** = 'RMP\_RESET'

Reset a selected ramp program

**RAMP\_ITERATIONS** = 'RMP\_OUT\_02\_'

Define how often a ramp program should be iterated

**RAMP\_PAUSE** = 'RMP\_PAUSE'

Pause a selected ramp program

**RAMP\_SELECT** = 'RMP\_SELECT\_'

Select a ramp program (target for all further ramp commands)

```

RAMP_SET = 'RMP_OUT_00_'
    Define parameters of a selected ramp program

RAMP_START = 'RMP_START'
    Start a selected ramp program

RAMP_STOP = 'RMP_STOP'
    Stop a running ramp program

START = 'START'
    Start temp control (pump and heating/cooling)

STOP = 'STOP'
    Stop temp control (pump and heating/cooling)

TEMP_SET_POINT = 'OUT_SP_00_'
    Define temperature set point

UPPER_TEMP = 'OUT_SP_04_'
    Define upper temp limit

build_str(param: str = '', terminator: str = '\n')
    Build a command string for sending to the device

```

#### Parameters

- **param** – Command's parameter given as string
- **terminator** – Command's terminator

**Returns** Command's string with a parameter and terminator

#### exception **LaudaProRp245eCommandError**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

Exception raised when an error is returned upon a command.

```

class LaudaProRp245eConfig(temp_set_point_init: Union[int, float] = 20.0, pump_init: int = 6, upper_temp:
    Union[int, float] = 80.0, lower_temp: Union[int, float] = - 55.0, com_time_out:
    Union[int, float] = 0, max_pump_level: int = 8, max_pr_number: int = 5,
    operation_mode: Union[int,
    hvl\_ccb.dev.lauda.LaudaProRp245eConfig.OperationModeEnum] =
    OperationModeEnum.AUTO, control_mode: Union[int,
    hvl\_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] =
    ExtControlModeEnum.INTERNAL)

```

Bases: object

Configuration for the Lauda RP245E circulation chiller.

```

class ExtControlModeEnum(value)

```

Bases: [enum.IntEnum](#)

Source for definition of external, controlled temperature (option 2, 3 and 4 are not available with current configuration of the Lauda RP245E, add-on hardware would required)

```

ANALOG = 2

```

```

ETH = 5

```

```

EXPT100 = 1

```

```

INTERNAL = 0

```

```

SERIAL = 3

```

```
USB = 4
```

**class OperationModeEnum**(*value*)  
Bases: `enum.IntEnum`

Operation Mode (Cooling OFF/Cooling On/AUTO - set to AUTO)

```
AUTO = 2
```

Automatically select heating/cooling

```
COOLOFF = 0
```

```
COOLON = 1
```

**clean\_values()** → *None*

```
com_time_out: Union[int, float] = 0
```

Communication time out (0 = OFF)

```
control_mode: Union[int, hv1_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = 0
```

**force\_value**(*fieldname*, *value*)  
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

```
is_configdataclass = True
```

**classmethod keys()** → `Sequence[str]`  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

```
lower_temp: Union[int, float] = -55.0
```

Lower temperature limit (safe for Galden HT135 cooling liquid)

```
max_pr_number: int = 5
```

Maximum number of ramp programs that can be stored in the memory of the chiller

```
max_pump_level: int = 8
```

Highest pump level of the chiller

```
operation_mode: Union[int, hv1_ccb.dev.lauda.LaudaProRp245eConfig.OperationModeEnum] = 2
```

**classmethod optional\_defaults()** → `Dict[str, object]`  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```
pump_init: int = 6
```

Default pump Level

**classmethod required\_keys()** → `Sequence[str]`  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**temp\_set\_point\_init:** `Union[int, float] = 20.0`

Default temperature set point

**upper\_temp:** `Union[int, float] = 80.0`

Upper temperature limit (safe for Galden HT135 cooling liquid)

**class** `LaudaProRp245eTcpCommunication(configuration)`

Bases: `hvl_ccb.comm.tcp.Tcp`

Implements the Communication Protocol for Lauda PRO RP245E TCP connection.

**close()** → *None*

Close the Lauda PRO RP245E TCP connection.

**static config\_cls()** → `Type[hvl_ccb.dev.lauda.LaudaProRp245eTcpCommunicationConfig]`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**open()** → *None*

Open the Lauda PRO RP245E TCP connection.

**Raises** `LaudaProRp245eCommandError` – if the connection fails.

**query\_command**(*command: hvl\_ccb.dev.lauda.LaudaProRp245eCommand, param: str = ""*) → *str*

Send and receive function. E.g. to be used when setting/changing device setting. :param *command*: first part of command string, defined in `LaudaProRp245eCommand` :param *param*: second part of command string, parameter (by default '') :return: *None*

**read()** → *str*

Receive value function. :return: reply from device as a string, the terminator, as well as the 'OK' stripped from the reply to make it directly useful as a value (e.g. in case the internal bath temperature is requested)

**write\_command**(*command: hvl\_ccb.dev.lauda.LaudaProRp245eCommand, param: str = ""*) → *None*

Send command function. :param *command*: first part of command string, defined in `LaudaProRp245eCommand` :param *param*: second part of command string, parameter (by default '') :return: *None*

**class** `LaudaProRp245eTcpCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], port: int = 54321, bufsize: int = 1024, wait_sec_pre_read_or_write: Union[int, float] = 0.005, terminator: str = '\n')`

Bases: `hvl_ccb.comm.tcp.TcpCommunicationConfig`

Configuration dataclass for `LaudaProRp245eTcpCommunication`.

**clean\_values()** → *None*

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod** `keys()` → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


**terminator:** str = '\r\n'

The terminator character

**wait\_sec\_pre\_read\_or\_write:** Union[int, float] = 0.005

Delay time between commands in seconds

### hvl\_ccb.dev.mbw973

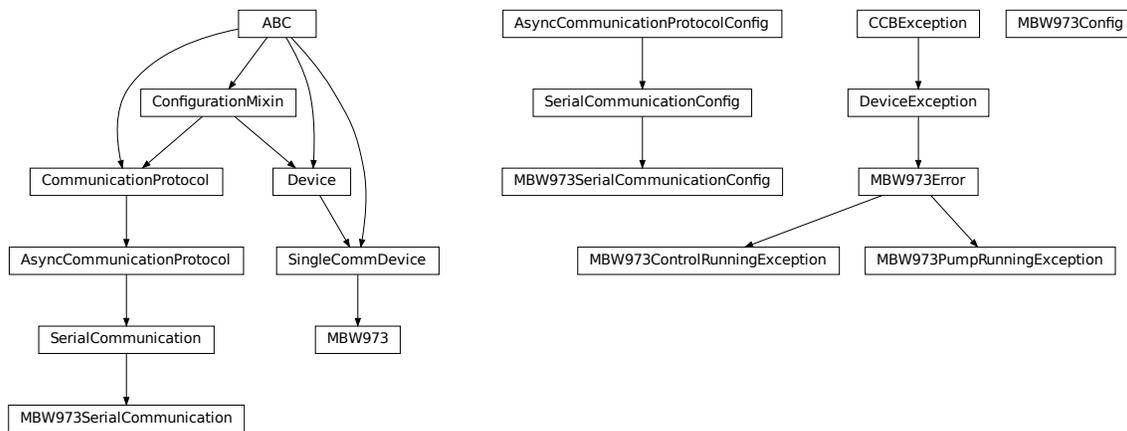

Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

**class MBW973**(com, dev\_config=None)

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

MBW 973 dew point mirror device class.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**is\_done()** → bool

Poll status of the dew point mirror and return True, if all measurements are done.

**Returns** True, if all measurements are done; False otherwise.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**read**(*cast\_type: typing.Type = <class 'str'>*)

Read value from *self.com* and cast to *cast\_type*. Raises *ValueError* if read text (*str*) is not convertible to *cast\_type*, e.g. to *float* or to *int*.

**Returns** Read value of *cast\_type* type.

**read\_float()** → float

Convenience wrapper for *self.read()*, with typing hint for return value.

**Returns** Read *float* value.

**read\_int()** → int

Convenience wrapper for *self.read()*, with typing hint for return value.

**Returns** Read *int* value.

**read\_measurements()** → Dict[str, float]

Read out measurement values and return them as a dictionary.

**Returns** Dictionary with values.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_measuring\_options**(*humidity: bool = True, sf6\_purity: bool = False*) → *None*

Send measuring options to the dew point mirror.

**Parameters**

- **humidity** – Perform humidity test or not?
- **sf6\_purity** – Perform SF6 purity test or not?

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**start()** → *None*

Start this device. Opens the communication protocol and retrieves the set measurement options from the device.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened.

**start\_control()** → *None*

Start dew point control to acquire a new value set.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**stop()** → *None*

Stop the device. Closes also the communication protocol.

**write**(*value*) → *None*

Send *value* to *self.com*.

**Parameters** **value** – Value to send, converted to *str*.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**class** **MBW973Config**(*polling\_interval: Union[int, float] = 2*)

Bases: object

Device configuration dataclass for MBW973.

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_interval: Union[int, float] = 2**

Polling period for *is\_done* status queries [in seconds].

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception MBW973ControlRunningException**

Bases: [hvl\\_ccb.dev.mbw973.MBW973Error](#)

Error indicating there is still a measurement running, and a new one cannot be started.

**exception MBW973Error**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

General error with the MBW973 dew point mirror device.

**exception MBW973PumpRunningException**

Bases: [hvl\\_ccb.dev.mbw973.MBW973Error](#)

Error indicating the pump of the dew point mirror is still recovering gas, unable to start a new measurement.

**class MBW973SerialCommunication**(*configuration*)

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for the MBW973 dew point mirror. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class MBW973SerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                     encoding_error_handling: str = 'strict',
                                     wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                     default_n_attempts_read_text_nonempty: int = 10, port: Union[str,
                                     NoneType] = None, baudrate: int = 9600, parity: Union[str,
                                     hvl_ccb.comm.serial.SerialCommunicationParity] =
                                     <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
                                     hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                     <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
                                     hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                     <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                     Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate:** `int = 9600`

Baudrate for MBW973 is 9600 baud

**bytesize:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8`

One byte is eight bits long

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** `Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'`

MBW973 does not use parity

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1`

MBW973 does use one stop bit


**terminator:** `bytes = b'\r'`

The terminator is only CR

**timeout:** `Union[int, float] = 3`

use 3 seconds timeout as default

## hvl\_ccb.dev.newport

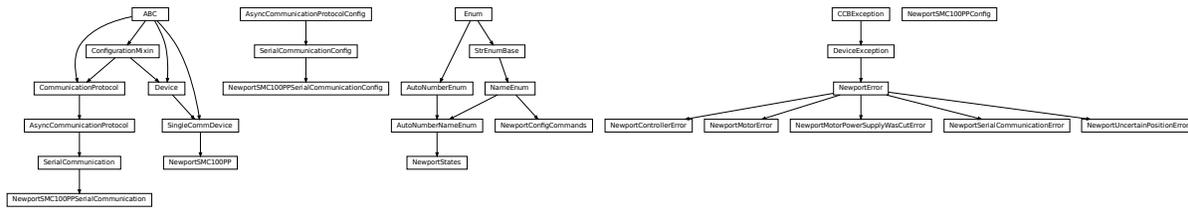

Device class for Newport SMC100PP stepper motor controller with serial communication.

The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/ff/smc100-single-axis-dc-or-stepper-motion-controller>

**class NewportConfigCommands** (*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl\_ccb.utils.enum.NameEnum*

Commands predefined by the communication protocol of the SMC100PP

**AC** = 'acceleration'

**BA** = 'backlash\_compensation'

**BH** = 'hysteresis\_compensation'

**FRM** = 'micro\_step\_per\_full\_step\_factor'

**FRS** = 'motion\_distance\_per\_full\_step'

**HT** = 'home\_search\_type'

**JR** = 'jerk\_time'

**OH** = 'home\_search\_velocity'

**OT** = 'home\_search\_timeout'

**QIL** = 'peak\_output\_current\_limit'

**SA** = 'rs485\_address'

**SL** = 'negative\_software\_limit'

**SR** = 'positive\_software\_limit'

**VA** = 'velocity'

**VB** = 'base\_velocity'

**ZX** = 'stage\_configuration'

**exception NewportControllerError**

Bases: *hvl\_ccb.dev.newport.NewportError*

Error with the Newport controller.

**exception NewportError**

Bases: *hvl\_ccb.dev.base.DeviceException*

General Exception for Newport Device

**exception NewportMotorError**Bases: *hvl\_ccb.dev.newport.NewportError*

Error with the Newport motor.

**exception NewportMotorPowerSupplyWasCutError**Bases: *hvl\_ccb.dev.newport.NewportError*

Error with the Newport motor after the power supply was cut and then restored, without interrupting the communication with the controller.

**class NewportSMC100PP**(*com, dev\_config=None*)Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Device class of the Newport motor controller SMC100PP

**class MotorErrors**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)Bases: *aenum.Enum*Possible motor errors reported by the motor during `get_state()`.**DC\_VOLTAGE\_TOO\_LOW** = 3**FOLLOWING\_ERROR** = 6**HOMING\_TIMEOUT** = 5**NED\_END\_OF\_TURN** = 11**OUTPUT\_POWER\_EXCEEDED** = 2**PEAK\_CURRENT\_LIMIT** = 9**POS\_END\_OF\_TURN** = 10**RMS\_CURRENT\_LIMIT** = 8**SHORT\_CIRCUIT** = 7**WRONG\_ESP\_STAGE** = 4**class StateMessages**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)Bases: *aenum.Enum*Possible messages returned by the controller on `get_state()` query.**CONFIG** = '14'**DISABLE\_FROM\_JOGGING** = '3E'**DISABLE\_FROM\_MOVING** = '3D'**DISABLE\_FROM\_READY** = '3C'**HOMING\_FROM\_RS232** = '1E'**HOMING\_FROM\_SMC** = '1F'**JOGGING\_FROM\_DISABLE** = '47'**JOGGING\_FROM\_READY** = '46'**MOVING** = '28'**NO\_REF\_ESP\_STAGE\_ERROR** = '10'**NO\_REF\_FROM\_CONFIG** = '0C'

```
NO_REF_FROM_DISABLED = '0D'  
NO_REF_FROM_HOMING = '0B'  
NO_REF_FROM_JOGGING = '11'  
NO_REF_FROM_MOVING = '0F'  
NO_REF_FROM_READY = '0E'  
NO_REF_FROM_RESET = '0A'  
READY_FROM_DISABLE = '34'  
READY_FROM_HOMING = '32'  
READY_FROM_JOGGING = '35'  
READY_FROM_MOVING = '33'
```

**States**

alias of `hvl_ccb.dev.newport.NewportStates`

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**exit\_configuration**(*add: Optional[int] = None*) → *None*

Exit the CONFIGURATION state and go back to the NOT REFERENCED state. All configuration parameters are saved to the device's memory.

**Parameters** *add* – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**get\_acceleration**(*add: Optional[int] = None*) → Union[int, float]

Leave the configuration state. The configuration parameters are saved to the device's memory.

**Parameters** *add* – controller address (1 to 31)

**Returns** acceleration (preset units/s<sup>2</sup>), value between 1e-6 and 1e12

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**get\_controller\_information**(*add: Optional[int] = None*) → str

Get information on the controller name and driver version

**Parameters** *add* – controller address (1 to 31)

**Returns** controller information

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_motor\_configuration**(*add: Optional[int] = None*) → Dict[str, float]

Query the motor configuration and returns it in a dictionary.

**Parameters** **add** – controller address (1 to 31)

**Returns** dictionary containing the motor's configuration

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_move\_duration**(*dist: Union[int, float], add: Optional[int] = None*) → float

Estimate the time necessary to move the motor of the specified distance.

**Parameters**

- **dist** – distance to travel
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_negative\_software\_limit**(*add: Optional[int] = None*) → Union[int, float]

Get the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

**Parameters** **add** – controller address (1 to 31)

**Returns** negative software limit (preset units), value between -1e12 and 0

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_position**(*add: Optional[int] = None*) → float

Returns the value of the current position.

**Parameters** **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportUncertainPositionError* – if the position is ambiguous

**get\_positive\_software\_limit**(*add: Optional[int] = None*) → Union[int, float]

Get the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

**Parameters** *add* – controller address (1 to 31)

**Returns** positive software limit (preset units), value between 0 and 1e12

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_state**(*add: int = None*) → *StateMessages*

Check on the motor errors and the controller state

**Parameters** *add* – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportMotorError* – if the motor reports an error

**Returns** state message from the device (member of *StateMessages*)

**go\_home**(*add: Optional[int] = None*) → *None*

Move the motor to its home position.

**Parameters** *add* – controller address (1 to 31), defaults to *self.address*

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**go\_to\_configuration**(*add: Optional[int] = None*) → *None*

This method is executed during *start()*. It can also be executed after a *reset()*. The controller is put in CONFIG state, where configuration parameters can be changed.

**Parameters** *add* – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**initialize**(*add: Optional[int] = None*) → *None*

Puts the controller from the NOT\_REF state to the READY state. Sends the motor to its “home” position.

**Parameters** *add* – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed

- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**move\_to\_absolute\_position**(*pos*: Union[int, float], *add*: Optional[int] = None) → None

Move the motor to the specified position.

#### Parameters

- **pos** – target absolute position (affected by the configured offset)
- **add** – controller address (1 to 31), defaults to self.address

#### Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**move\_to\_relative\_position**(*pos*: Union[int, float], *add*: Optional[int] = None) → None

Move the motor of the specified distance.

#### Parameters

- **pos** – distance to travel (the sign gives the direction)
- **add** – controller address (1 to 31), defaults to self.address

#### Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**reset**(*add*: Optional[int] = None) → None

Resets the controller, equivalent to a power-up. This puts the controller back to NOT REFERENCED state, which is necessary for configuring the controller.

**Parameters** **add** – controller address (1 to 31)

#### Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**set\_acceleration**(*acc*: Union[int, float], *add*: Optional[int] = None) → None

Leave the configuration state. The configuration parameters are saved to the device's memory.

#### Parameters

- **acc** – acceleration (preset units/s<sup>2</sup>), value between 1e-6 and 1e12
- **add** – controller address (1 to 31)

#### Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**set\_motor\_configuration**(*add: Optional[int] = None, config: Optional[dict] = None*) → *None*

Set the motor configuration. The motor must be in CONFIG state.

**Parameters**

- **add** – controller address (1 to 31)
- **config** – dictionary containing the motor's configuration

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**set\_negative\_software\_limit**(*lim: Union[int, float], add: Optional[int] = None*) → *None*

Set the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

**Parameters**

- **lim** – negative software limit (preset units), value between -1e12 and 0
- **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**set\_positive\_software\_limit**(*lim: Union[int, float], add: Optional[int] = None*) → *None*

Set the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

**Parameters**

- **lim** – positive software limit (preset units), value between 0 and 1e12
- **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**start()**

Opens the communication protocol and applies the config.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened

**stop()** → *None*

Stop the device. Close the communication protocol.

**stop\_motion**(*add: Optional[int] = None*) → *None*

Stop a move in progress by decelerating the positioner immediately with the configured acceleration until it stops. If a controller address is provided, stops a move in progress on this controller, else stops the moves on all controllers.

**Parameters** **add** – controller address (1 to 31)

**Raises**

- **`SerialCommunicationIOError`** – if the com is closed
- **`NewportSerialCommunicationError`** – if an unexpected answer is obtained
- **`NewportControllerError`** – if the controller reports an error

**`wait_until_motor_initialized`**(*add: Optional[int] = None*) → *None*

Wait until the motor leaves the HOMING state (at which point it should have arrived to the home position).

**Parameters** *add* – controller address (1 to 31)

**Raises**

- **`SerialCommunicationIOError`** – if the com is closed
- **`NewportSerialCommunicationError`** – if an unexpected answer is obtained
- **`NewportControllerError`** – if the controller reports an error

```
class NewportSMC100PPConfig(address: int = 1, user_position_offset: Union[int, float] = 23.987,
                           screw_scaling: Union[int, float] = 1, exit_configuration_wait_sec: Union[int,
                           float] = 5, move_wait_sec: Union[int, float] = 1, acceleration: Union[int, float]
                           = 10, backlash_compensation: Union[int, float] = 0, hysteresis_compensation:
                           Union[int, float] = 0.015, micro_step_per_full_step_factor: int = 100,
                           motion_distance_per_full_step: Union[int, float] = 0.01, home_search_type:
                           Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] =
                           HomeSearch.HomeSwitch, jerk_time: Union[int, float] = 0.04,
                           home_search_velocity: Union[int, float] = 4, home_search_timeout: Union[int,
                           float] = 27.5, home_search_polling_interval: Union[int, float] = 1,
                           peak_output_current_limit: Union[int, float] = 0.4, rs485_address: int = 2,
                           negative_software_limit: Union[int, float] = - 23.5, positive_software_limit:
                           Union[int, float] = 25, velocity: Union[int, float] = 4, base_velocity: Union[int,
                           float] = 0, stage_configuration: Union[int,
                           hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageConfig] =
                           EspStageConfig.EnableEspStageCheck)
```

Bases: object

Configuration dataclass for the Newport motor controller SMC100PP.

```
class EspStageConfig(value=<no_arg>, names=None, module=None, type=None, start=1,
                    boundary=None)
```

Bases: aenum.IntEnum

Different configurations to check or not the motor configuration upon power-up.

**DisableEspStageCheck** = 1

**EnableEspStageCheck** = 3

**UpdateEspStageInfo** = 2

```
class HomeSearch(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: aenum.IntEnum

Different methods for the motor to search its home position during initialization.

**CurrentPosition** = 1

**EndOfRunSwitch** = 4

**EndOfRunSwitch\_and\_Index** = 3

**HomeSwitch** = 2

```
    HomeSwitch_and_Index = 0
acceleration: Union[int, float] = 10
address: int = 1
backlash_compensation: Union[int, float] = 0
base_velocity: Union[int, float] = 0
clean_values()
exit_configuration_wait_sec: Union[int, float] = 5
force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing
    after value has been forced on fieldname.
```

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

```
home_search_polling_interval: Union[int, float] = 1
home_search_timeout: Union[int, float] = 27.5
home_search_type: Union[int, hvl\_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch]
= 2
home_search_velocity: Union[int, float] = 4
hysteresis_compensation: Union[int, float] = 0.015
is_configdataclass = True
jerk_time: Union[int, float] = 0.04
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.
micro_step_per_full_step_factor: int = 100
motion_distance_per_full_step: Union[int, float] = 0.01
property motor_config: Dict[str, float]
    Gather the configuration parameters of the motor into a dictionary.

    Returns dict containing the configuration parameters of the motor
move_wait_sec: Union[int, float] = 1
negative_software_limit: Union[int, float] = -23.5
classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
    on instantiation.

    Returns a list of strings containing all optional keys.
peak_output_current_limit: Union[int, float] = 0.4
positive_software_limit: Union[int, float] = 25
```

`post_force_value`(*fieldname*, *value*)

**classmethod** `required_keys`() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

`rs485_address`: int = 2

`screw_scaling`: Union[int, float] = 1

`stage_configuration`: Union[int, [hvl\\_ccb.dev.newport NewportSMC100PPConfig.EspStageConfig](#)] = 3

`user_position_offset`: Union[int, float] = 23.987

`velocity`: Union[int, float] = 4

**class** `NewportSMC100PPSerialCommunication`(*configuration*)

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for NewportSMC100 controller. Already predefines device-specific protocol parameters in config.

**class** `ControllerErrors`(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `aenum.Enum`

Possible controller errors with values as returned by the device in response to sent commands.

`ADDR_INCORRECT` = 'B'

`CMD_EXEC_ERROR` = 'V'

`CMD_NOT_ALLOWED` = 'D'

`CMD_NOT_ALLOWED_CC` = 'X'

`CMD_NOT_ALLOWED_CONFIGURATION` = 'I'

`CMD_NOT_ALLOWED_DISABLE` = 'J'

`CMD_NOT_ALLOWED_HOMING` = 'L'

`CMD_NOT_ALLOWED_MOVING` = 'M'

`CMD_NOT_ALLOWED_NOT_REFERENCED` = 'H'

`CMD_NOT_ALLOWED_PP` = 'W'

`CMD_NOT_ALLOWED_READY` = 'K'

`CODE_OR_ADDR_INVALID` = 'A'

`COM_TIMEOUT` = 'S'

`DISPLACEMENT_OUT_OF_LIMIT` = 'G'

`EEPROM_ACCESS_ERROR` = 'U'

`ESP_STAGE_NAME_INVALID` = 'F'

`HOME_STARTED` = 'E'

`NO_ERROR` = '@'

`PARAM_MISSING_OR_INVALID` = 'C'

`POSITION_OUT_OF_LIMIT = 'N'`

`check_for_error(add: int) → None`

Ask the Newport controller for the last error it recorded.

This method is called after every command or query.

**Parameters** `add` – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

`static config_cls()`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

`query(add: int, cmd: str, param: Optional[Union[int, float, str]] = None) → str`

Send a query to the controller, read the answer, and check for errors. The prefix `add+cmd` is removed from the answer.

**Parameters**

- `add` – the controller address (1 to 31)
- `cmd` – the command to be sent
- `param` – optional parameter (int/float/str) appended to the command

**Returns** the answer from the device without the prefix

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

`query_multiple(add: int, cmd: str, prefixes: List[str]) → List[str]`

Send a query to the controller, read the answers, and check for errors. The prefixes are removed from the answers.

**Parameters**

- `add` – the controller address (1 to 31)
- `cmd` – the command to be sent
- `prefixes` – prefixes of each line expected in the answer

**Returns** list of answers from the device without prefix

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**read\_text()** → str

Read one line of text from the serial port, and check for presence of a null char which indicates that the motor power supply was cut and then restored. The input buffer may hold additional data afterwards, since only one line is read.

This method uses *self.access\_lock* to ensure thread-safety.

**Returns** String read from the serial port; '' if there was nothing to read.

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **NewportMotorPowerSupplyWasCutError** – if a null char is read

**send\_command**(*add: int, cmd: str, param: Optional[Union[int, float, str]] = None*) → None

Send a command to the controller, and check for errors.

**Parameters**

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**send\_stop**(*add: int*) → None

Send the general stop ST command to the controller, and check for errors.

**Parameters** **add** – the controller address (1 to 31)

**Returns** ControllerErrors reported by Newport Controller

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained

```
class NewportSMC100PPSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'ascii',
                                                encoding_error_handling: str = 'replace',
                                                wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                                default_n_attempts_read_text_nonempty: int = 10,
                                                port: Union[str, NoneType] = None, baudrate: int =
                                                57600, parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 10)
```

Bases: *hvl\_ccb.comm.serial.SerialCommunicationConfig*

**baudrate:** `int = 57600`

Baudrate for NewportSMC100 controller is 57600 baud

**bytesize:** `Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8`

NewportSMC100 controller uses 8 bits for one data byte

**encoding:** `str = 'ascii'`

use ASCII as de-/encoding, cf. the manual

**encoding\_error\_handling:** `str = 'replace'`

replace bytes with `instead` of raising utf-8 exception when decoding fails

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod** `keys()` → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod** `optional_defaults()` → `Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** `Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'`

NewportSMC100 controller does not use parity

**classmethod** `required_keys()` → `Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** `Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1`

NewportSMC100 controller uses one stop bit

**terminator:** `bytes = b'\r\n'`

The terminator is CR/LF

**timeout:** `Union[int, float] = 10`

use 10 seconds timeout as default

**exception** `NewportSerialCommunicationError`

Bases: [hvl\\_ccb.dev.newport.NewportError](#)

Communication error with the Newport controller.

**class** `NewportStates`(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [hvl\\_ccb.utils.enum.AutoNumberNameEnum](#)

States of the Newport controller. Certain commands are allowed only in certain states.

**CONFIG** = 3

**DISABLE** = 6


**HOMING** = 2  
**JOGGING** = 7  
**MOVING** = 5  
**NO\_REF** = 1  
**READY** = 4

**exception NewportUncertainPositionError**

Bases: *hvl\_ccb.dev.newport.NewportError*

Error with the position of the Newport motor.

**hvl\_ccb.dev.pfeiffer\_tpg**

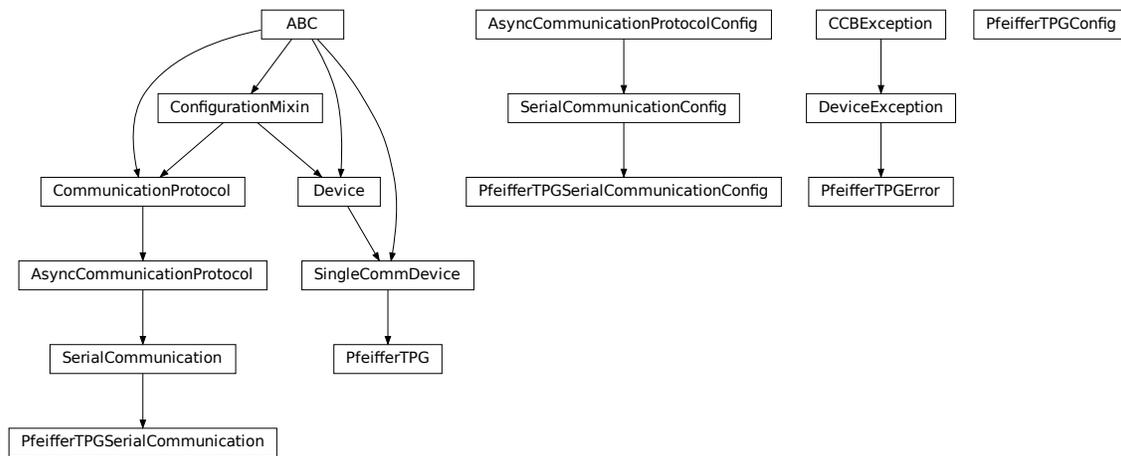

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/measurement/activeline/controllers/>

**class PfeifferTPG**(*com, dev\_config=None*)

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Pfeiffer TPG control unit device class

**class PressureUnits**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl\_ccb.utils.enum.NameEnum*

Enum of available pressure units for the digital display. “0” corresponds either to bar or to mbar depending on the TPG model. In case of doubt, the unit is visible on the digital display.

**Micron** = 3  
**Pascal** = 2  
**Torr** = 1

```
Volt = 5
bar = 0
hPascal = 4
mbar = 0
class SensorStatus(value)
    Bases: enum.IntEnum
    An enumeration.
    Identification_error = 6
    No_sensor = 5
    Ok = 0
    Overrange = 2
    Sensor_error = 3
    Sensor_off = 4
    Underrange = 1
class SensorTypes(value)
    Bases: enum.Enum
    An enumeration.
    CMR = 4
    IKR = 2
    IKR11 = 2
    IKR9 = 2
    IMR = 5
    None = 7
    PBR = 6
    PKR = 3
    TPR = 1
    noSENSOR = 7
    noSen = 7
static config_cls()
    Return the default configdataclass class.
    Returns a reference to the default configdataclass class
static default_com_cls()
    Get the class for the default communication protocol used with this device.
    Returns the type of the standard communication protocol for this device
get_full_scale_mbar() → List[Union[int, float]]
    Get the full scale range of the attached sensors
    Returns full scale range values in mbar, like [0.01, 1, 0.1, 1000, 50000, 10]
    Raises
```

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**get\_full\_scale\_unitless()** → List[int]

Get the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

**Returns** list of full scale range values, like *[0, 1, 3, 3, 2, 0]*

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**identify\_sensors()** → *None*

Send identification request TID to sensors on all channels.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**measure(channel: int)** → Tuple[str, float]

Get the status and measurement of one sensor

**Parameters** **channel** – int channel on which the sensor is connected, with  $1 \leq \text{channel} \leq \text{number\_of\_sensors}$

**Returns** measured value as float if measurement successful, sensor status as string if not

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**measure\_all()** → List[Tuple[str, float]]

Get the status and measurement of all sensors (this command is not available on all models)

**Returns** list of measured values as float if measurements successful, and or sensor status as strings if not

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**property number\_of\_sensors**

**set\_display\_unit(unit: Union[str, hvl\_ccb.dev.pfeiffer\_tpg.PfeifferTPG.PressureUnits])** → *None*

Set the unit in which the measurements are shown on the display.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**set\_full\_scale\_mbar(fsr: List[Union[int, float]])** → *None*

Set the full scale range of the attached sensors (in unit mbar)

**Parameters** **fsr** – full scale range values in mbar, for example *[0.01, 1000]*

**Raises**

- **`SerialCommunicationIOError`** – when communication port is not opened
- **`PfeifferTPGError`** – if command fails

**set\_full\_scale\_unitless**(*fsr*: List[int]) → None

Set the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

**Parameters** *fsr* – list of full scale range values, like [0, 1, 3, 3, 2, 0]

**Raises**

- **`SerialCommunicationIOError`** – when communication port is not opened
- **`PfeifferTPGError`** – if command fails

**start**() → None

Start this device. Opens the communication protocol, and identify the sensors.

**Raises** **`SerialCommunicationIOError`** – when communication port cannot be opened

**stop**() → None

Stop the device. Closes also the communication protocol.

**property unit**

The pressure unit of readings is always mbar, regardless of the display unit.

**class PfeifferTPGConfig**(*model*: Union[str, hvl\_ccb.dev.pfeiffer\_tpg.PfeifferTPGConfig.Model] = Model.TPG25xA)

Bases: object

Device configuration dataclass for Pfeiffer TPG controllers.

**class Model**(*value*=<no\_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: `hvl_ccb.utils.enum.NameEnum`

An enumeration.

**TPG25xA** = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

**TPGx6x** = {0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8, 50000: 9}

**is\_valid\_scale\_range\_reversed\_str**(*v*: str) → bool

Check if given string represents a valid reversed scale range of a model.

**Parameters** *v* – Reversed scale range string.

**Returns** *True* if valid, *False* otherwise.

**clean\_values**()

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass** = True

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**model:** Union[str, [hvl\\_ccb.dev.pfeiffer\\_tpg.PfeifferTPGConfig.Model](#)] = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception PfeifferTPGError**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

Error with the Pfeiffer TPG Controller.

**class PfeifferTPGSerialCommunication(configuration)**

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for Pfeiffer TPG controllers. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query(cmd: str)** → str

Send a query, then read and returns the first line from the com port.

**Parameters cmd** – query message to send to the device

**Returns** first line read on the com

**Raises**

- [SerialCommunicationIOError](#) – when communication port is not opened
- [PfeifferTPGError](#) – if the device does not acknowledge the command or if the answer from the device is empty

**send\_command(cmd: str)** → None

Send a command to the device and check for acknowledgement.

**Parameters cmd** – command to send to the device

**Raises**

- [SerialCommunicationIOError](#) – when communication port is not opened
- [PfeifferTPGError](#) – if the answer from the device differs from the expected acknowledgement character 'chr(6)'.

```
class PfeifferTPGSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
    encoding_error_handling: str = 'strict',
    wait_sec_read_text_nonempty: Union[int, float] = 0.5,
    default_n_attempts_read_text_nonempty: int = 10, port:
    Union[str, NoneType] = None, baudrate: int = 9600, parity:
    Union[str,
    hvl_ccb.comm.serial.SerialCommunicationParity] =
    <SerialCommunicationParity.NONE: 'N'>, stopbits:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationStopbits] =
    <SerialCommunicationStopbits.ONE: 1>, bytesize:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationBytesize] =
    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
    Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate:** `int = 9600`

Baudrate for Pfeiffer TPG controllers is 9600 baud

**bytesize:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8`

One byte is eight bits long

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** `Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'`

Pfeiffer TPG controllers do not use parity

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1`

Pfeiffer TPG controllers use one stop bit

**terminator:** `bytes = b'\r\n'`

The terminator is <CR><LF>

```
timeout: Union[int, float] = 3
    use 3 seconds timeout as default
```

### hvl\_ccb.dev.picotech\_pt104


**NOTE:** PicoSDK Python wrappers already on import attempt to load the PicoSDK library; thus, the API docs can only be generated in a system with the latter installed and are by default disabled.

To build the API docs for this submodule locally edit the docs/hvl\_ccb.dev.picotech\_pt104.rst file to remove the .. code-block:: directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.picotech_pt104
    :parts: 1

.. automodule:: hvl_ccb.dev.picotech_pt104
    :members:
    :undoc-members:
    :show-inheritance:
```

### hvl\_ccb.dev.rs\_rto1024

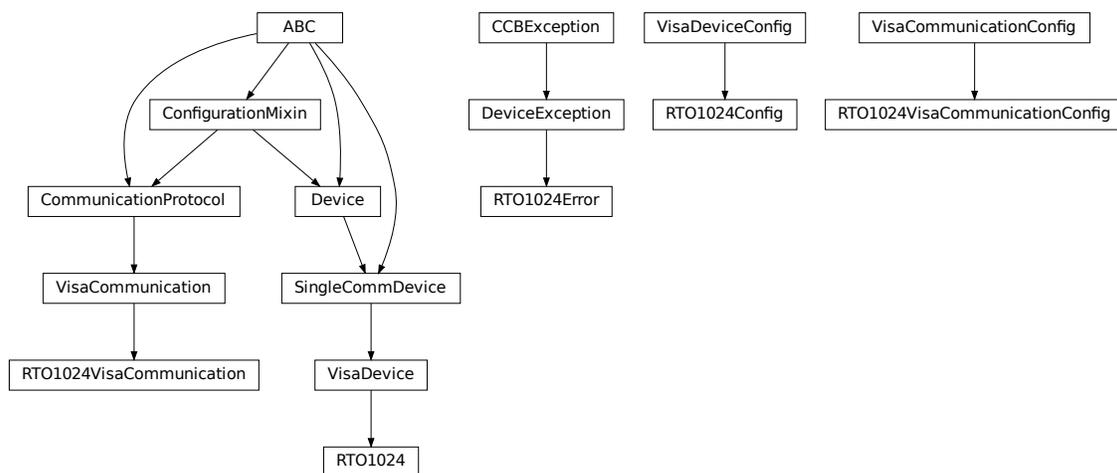

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

```
class RTO1024(com: Union[hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunication,
    hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunicationConfig, dict], dev_config:
    Union[hvl_ccb.dev.rs_rto1024.RTO1024Config, dict])
```

Bases: *hvl\_ccb.dev.visa.VisaDevice*

Device class for the Rhode & Schwarz RTO 1024 oscilloscope.

```
class TriggerModes(value=<no_arg>, names=None, module=None, type=None, start=1,
    boundary=None)
```

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

Enumeration for the three available trigger modes.

**AUTO = 1**

**FREERUN = 3**

**NORMAL = 2**

**classmethod names()**

Returns a list of the available trigger modes. :return: list of strings

**activate\_measurements**(*meas\_n: int, source: str, measurements: List[str], category: str = 'AMPTime'*)

Activate the list of 'measurements' of the waveform 'source' in the measurement box number 'meas\_n'. The list 'measurements' starts with the main measurement and continues with additional measurements of the same 'category'.

**Parameters**

- **meas\_n** – measurement number 1..8
- **source** – measurement source, for example C1W1
- **measurements** – list of measurements, the first one will be the main measurement.
- **category** – the category of measurements, by default AMPTime

**backup\_waveform**(*filename: str*) → *None*

Backup a waveform file from the standard directory specified in the device configuration to the standard backup destination specified in the device configuration. The filename has to be specified without .bin or path.

**Parameters filename** – The waveform filename without extension and path

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns** the VisaCommunication class

**file\_copy**(*source: str, destination: str*) → *None*

Copy a file from one destination to another on the oscilloscope drive. If the destination file already exists, it is overwritten without notice.

**Parameters**

- **source** – absolute path to the source file on the DSO filesystem
- **destination** – absolute path to the destination file on the DSO filesystem

**Raises** *RT01024Error* – if the operation did not complete

**get\_acquire\_length**() → float

Gets the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]

**Returns** the time for one acquisition. Range: 250e-12 ... 500 [s]

**get\_channel\_offset**(*channel: int*) → float

Gets the voltage offset of the indicated channel.

**Parameters** **channel** – is the channel number (1..4)

**Returns** channel offset voltage in V (value between -1 and 1)

**get\_channel\_position**(*channel: int*) → float

Gets the vertical position of the indicated channel.

**Parameters** **channel** – is the channel number (1..4)

**Returns** channel position in div (value between -5 and 5)

**get\_channel\_range**(*channel: int*) → float

Queries the channel range in V.

**Parameters** **channel** – is the input channel (1..4)

**Returns** channel range in V

**get\_channel\_scale**(*channel: int*) → float

Queries the channel scale in V/div.

**Parameters** **channel** – is the input channel (1..4)

**Returns** channel scale in V/div

**get\_channel\_state**(*channel: int*) → bool

Queries if the channel is active or not.

**Parameters** **channel** – is the input channel (1..4)

**Returns** True if active, else False

**get\_reference\_point**() → int

Gets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 ... 100 [%]
- Increment: 1 [%]

**Returns** the reference in %

**get\_repetitions**() → int

Get the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 ... 16777215
- Increment: 10
- \*RST = 1

**Returns** the number of waveforms to acquire

**get\_timestamps**() → List[float]

Gets the timestamps of all recorded frames in the history and returns them as a list of floats.

**Returns** list of timestamps in [s]

**Raises** *RT01024Error* – if the timestamps are invalid

**list\_directory**(*path: str*) → List[Tuple[str, str, int]]

List the contents of a given directory on the oscilloscope filesystem.

**Parameters** **path** – is the path to a folder

**Returns** a list of filenames in the given folder

**load\_configuration**(*filename: str*) → *None*

Load current settings from a configuration file. The filename has to be specified without base directory and '.dff' extension.

**Information from the manual** *ReCaLl* calls up the instrument settings from an intermediate memory identified by the specified number. The instrument settings can be stored to this memory using the command *\*SAV* with the associated number. It also activates the instrument settings which are stored in a file and loaded using *MMEMemory:LOAD:STATe* .

**Parameters filename** – is the name of the settings file without path and extension

**local\_display**(*state: bool*) → *None*

Enable or disable local display of the scope.

**Parameters state** – is the desired local display state

**prepare\_ultra\_segmentation**() → *None*

Make ready for a new acquisition in ultra segmentation mode. This function does one acquisition without ultra segmentation to clear the history and prepare for a new measurement.

**read\_measurement**(*meas\_n: int, name: str*) → float

**Parameters**

- **meas\_n** – measurement number 1..8
- **name** – measurement name, for example “MAX”

**Returns** measured value

**run\_continuous\_acquisition**() → *None*

Start acquiring continuously.

**run\_single\_acquisition**() → *None*

Start a single or Nx acquisition.

**save\_configuration**(*filename: str*) → *None*

Save the current oscilloscope settings to a file. The filename has to be specified without path and '.dff' extension, the file will be saved to the configured settings directory.

**Information from the manual** *SAVe* stores the current instrument settings under the specified number in an intermediate memory. The settings can be recalled using the command *\*RCL* with the associated number. To transfer the stored instrument settings to a file, use *MMEMemory:STORe:STATe* .

**Parameters filename** – is the name of the settings file without path and extension

**save\_waveform\_history**(*filename: str, channel: int, waveform: int = 1*) → *None*

Save the history of one channel and one waveform to a .bin file. This function is used after an acquisition using sequence trigger mode (with or without ultra segmentation) was performed.

**Parameters**

- **filename** – is the name (without extension) of the file
- **channel** – is the channel number
- **waveform** – is the waveform number (typically 1)

**Raises** *RT01024Error* – if storing waveform times out

**set\_acquire\_length**(*timerange: float*) → *None*

Defines the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]
- \*RST = 0.5 [s]

**Parameters timerange** – is the time for one acquisition. Range: 250e-12 ... 500 [s]

**set\_channel\_offset**(*channel: int, offset: float*) → *None*

Sets the voltage offset of the indicated channel.

- Range: Dependent on the channel scale and coupling [V]
- Increment: Minimum 0.001 [V], may be higher depending on the channel scale and coupling
- \*RST = 0

#### Parameters

- **channel** – is the channel number (1..4)
- **offset** – Offset voltage. Positive values move the waveform down, negative values move it up.

**set\_channel\_position**(*channel: int, position: float*) → *None*

Sets the vertical position of the indicated channel as a graphical value.

- Range: -5.0 ... 5.0 [div]
- Increment: 0.02
- \*RST = 0

#### Parameters

- **channel** – is the channel number (1..4)
- **position** – is the position. Positive values move the waveform up, negative values move it down.

**set\_channel\_range**(*channel: int, v\_range: float*) → *None*

Sets the voltage range across the 10 vertical divisions of the diagram. Use the command alternatively instead of set\_channel\_scale.

- Range for range: Depends on attenuation factors and coupling. With 1:1 probe and external attenuations and 50 input coupling, the range is 10 mV to 10 V. For 1 M input coupling, it is 10 mV to 100 V. If the probe and/or external attenuation is changed, multiply the range values by the attenuation factors.
- Increment: 0.01
- \*RST = 0.5

#### Parameters

- **channel** – is the channel number (1..4)
- **v\_range** – is the vertical range [V]

**set\_channel\_scale**(*channel: int, scale: float*) → *None*

Sets the vertical scale for the indicated channel. The scale value is given in volts per division.

- Range for scale: depends on attenuation factor and coupling. With 1:1 probe and external attenuations and 50 input coupling, the vertical scale (input sensitivity) is 1 mV/div to 1 V/div. For 1 M input coupling, it is 1 mV/div to 10 V/div. If the probe and/or external attenuation is changed, multiply the values by the attenuation factors to get the actual scale range.
- Increment: 1e-3
- \*RST = 0.05

See also: `set_channel_range`

**Parameters**

- **channel** – is the channel number (1..4)
- **scale** – is the vertical scaling [V/div]

`set_channel_state(channel: int, state: bool) → None`

Switches the channel signal on or off.

**Parameters**

- **channel** – is the input channel (1..4)
- **state** – is True for on, False for off

`set_reference_point(percentage: int) → None`

Sets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 ... 100 [%]
- Increment: 1 [%]
- \*RST = 50 [%]

**Parameters percentage** – is the reference in %

`set_repetitions(number: int) → None`

Set the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 ... 16777215
- Increment: 10
- \*RST = 1

**Parameters number** – is the number of waveforms to acquire

`set_trigger_level(channel: int, level: float, event_type: int = 1) → None`

Sets the trigger level for the specified event and source.

- Range: -10 to 10 V
- Increment: 1e-3 V
- \*RST = 0 V

**Parameters**

- **channel** – indicates the trigger source.
  - 1..4 = channel 1 to 4, available for all event types 1..3

- 5 = external trigger input on the rear panel for analog signals, available for A-event type = 1
- 6..9 = not available
- **level** – is the voltage for the trigger level in [V].
- **event\_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

**set\_trigger\_mode**(*mode: Union[str, hvl\_ccb.dev.rs\_rto1024.RTO1024.TriggerModes]*) → *None*  
Sets the trigger mode which determines the behavior of the instrument if no trigger occurs.

**Parameters mode** – is either auto, normal, or freerun.

**Raises *RT01024Error*** – if an invalid triggermode is selected

**set\_trigger\_source**(*channel: int, event\_type: int = 1*) → *None*  
Set the trigger (Event A) source channel.

**Parameters**

- **channel** – is the channel number (1..4)
- **event\_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

**start**() → *None*

Start the RTO1024 oscilloscope and bring it into a defined state and remote mode.

**stop**() → *None*

Stop the RTO1024 oscilloscope, reset events and close communication. Brings back the device to a state where local operation is possible.

**stop\_acquisition**() → *None*  
Stop any acquisition.

**class RT01024Config**(*waveforms\_path: str, settings\_path: str, backup\_path: str, spoll\_interval: Union[int, float] = 0.5, spoll\_start\_delay: Union[int, float] = 2, command\_timeout\_seconds: Union[int, float] = 60, wait\_sec\_short\_pause: Union[int, float] = 0.1, wait\_sec\_enable\_history: Union[int, float] = 1, wait\_sec\_post\_acquisition\_start: Union[int, float] = 2*)

Bases: *hvl\_ccb.dev.visa.VisaDeviceConfig*, *hvl\_ccb.dev.rs\_rto1024.\_RT01024ConfigDefaultsBase*, *hvl\_ccb.dev.rs\_rto1024.\_RT01024ConfigBase*

Configdataclass for the RTO1024 device.

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod** `required_keys()` → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception** `RT01024Error`

Bases: `hvl_ccb.dev.base.DeviceException`

**class** `RT01024VisaCommunication(configuration)`

Bases: `hvl_ccb.comm.visa.VisaCommunication`

Specialization of `VisaCommunication` for the RT01024 oscilloscope

**static** `config_cls()`

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class** `RT01024VisaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], interface_type: Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType] = InterfaceType.TCPIP_INSTR, board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = '')`

Bases: `hvl_ccb.comm.visa.VisaCommunicationConfig`

Configuration dataclass for `VisaCommunication` with specifications for the RT01024 device class.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**interface\_type: Union[str, hvl\_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]**  
**= 2**

Interface type of the VISA connection, being one of `InterfaceType`.

**classmethod** `keys()` → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod** `optional_defaults()` → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.


**Returns** a list of strings containing all optional keys.

**classmethod** `required_keys()` → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

## hvl\_ccb.dev.se\_ils2t

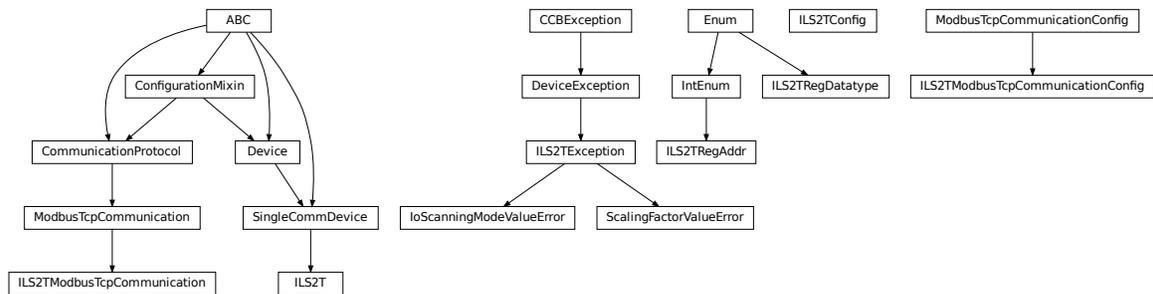

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

```
class ILS2T(com, dev_config=None)
```

Bases: `hvl_ccb.dev.base.SingleCommDevice`

Schneider Electric ILS2T stepper drive class.

```
ACTION_JOG_VALUE = 0
```

The single action value for `ILS2T.Mode.JOG`

```
class ActionsPtp(value)
```

Bases: `enum.IntEnum`

Allowed actions in the point to point mode (`ILS2T.Mode.PTP`).

```
ABSOLUTE_POSITION = 0
```

```
RELATIVE_POSITION_MOTOR = 2
```

```
RELATIVE_POSITION_TARGET = 1
```

```
DEFAULT_IO_SCANNING_CONTROL_VALUES = {'action': 2, 'continue_after_stop_cu': 0,
'disable_driver_di': 0, 'enable_driver_en': 0, 'execute_stop_sh': 0,
'fault_reset_fr': 0, 'mode': 3, 'quick_stop_qs': 0, 'ref_16': 1500, 'ref_32':
0, 'reset_stop_ch': 0}
```

Default IO Scanning control mode values

```
class Mode(value)
```

Bases: `enum.IntEnum`

ILS2T device modes

```
JOG = 1
```

```
PTP = 3
```

```
class Ref16Jog(value)
```

Bases: `enum.Flag`

Allowed values for ILS2T ref\_16 register (the shown values are the integer representation of the bits), all in Jog mode = 1

```
FAST = 4
```

```
NEG = 2
```

```
NEG_FAST = 6
```

```
NONE = 0
```

POS = 1

POS\_FAST = 5

**RegAddr**

Modbus Register Adresses

alias of *hvl\_ccb.dev.se\_ils2t.ILS2TRegAddr*

**RegDatatype**

Modbus Register Datatypes

alias of *hvl\_ccb.dev.se\_ils2t.ILS2TRegDatatype*

**class State(value)**

Bases: `enum.IntEnum`

State machine status values

ON = 6

QUICKSTOP = 7

READY = 4

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**disable(log\_warn: bool = True, wait\_sec\_max: Optional[int] = None) → bool**

Disable the driver of the stepper motor and enable the brake.

Note: the driver cannot be disabled if the motor is still running.

**Parameters**

- **log\_warn** – if log a warning in case the motor cannot be disabled.
- **wait\_sec\_max** – maximal wait time for the motor to stop running and to disable it; by default, with *None*, use a config value

**Returns** *True* if disable request could and was sent, *False* otherwise.

**do\_ioscanning\_write(\*\*kwargs: int) → None**

Perform a write operation using IO Scanning mode.

**Parameters** **kwargs** – Keyword-argument list with options to send, remaining are taken from the defaults.

**enable() → None**

Enable the driver of the stepper motor and disable the brake.

**execute\_absolute\_position(position: int) → bool**

Execute a absolute position change, i.e. enable motor, perform absolute position change, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

**Parameters** **position** – absolute position of motor in user defined steps.

**Returns** *True* if actual position is as expected, *False* otherwise.

**execute\_relative\_step**(*steps: int*) → bool

Execute a relative step, i.e. enable motor, perform relative steps, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

**Parameters** **steps** – Number of steps.

**Returns** *True* if actual position is as expected, *False* otherwise.

**get\_dc\_volt**() → float

Read the DC supply voltage of the motor.

**Returns** DC input voltage.

**get\_error\_code**() → Dict[int, Dict[str, Any]]

Read all messages in fault memory. Will read the full error message and return the decoded values. At the end the fault memory of the motor will be deleted. In addition, `reset_error` is called to re-enable the motor for operation.

**Returns** Dictionary with all information

**get\_position**() → int

Read the position of the drive and store into status.

**Returns** Position step value

**get\_status**() → Dict[str, int]

Perform an IO Scanning read and return the status of the motor.

**Returns** dict with status information.

**get\_temperature**() → int

Read the temperature of the motor.

**Returns** Temperature in degrees Celsius.

**jog\_run**(*direction: bool = True, fast: bool = False*) → *None*

Slowly turn the motor in positive direction.

**jog\_stop**() → *None*

Stop turning the motor in Jog mode.

**quickstop**() → *None*

Stops the motor with high deceleration rate and falls into error state. Reset with `reset_error` to recover into normal state.

**reset\_error**() → *None*

Resets the motor into normal state after quick stop or another error occurred.

**set\_jog\_speed**(*slow: int = 60, fast: int = 180*) → *None*

Set the speed for jog mode. Default values correspond to startup values of the motor.

**Parameters**

- **slow** – RPM for slow jog mode.
- **fast** – RPM for fast jog mode.

**set\_max\_acceleration**(*rpm\_minute: int*) → *None*

Set the maximum acceleration of the motor.

**Parameters** **rpm\_minute** – revolution per minute per minute

**set\_max\_deceleration**(*rpm\_minute: int*) → *None*

Set the maximum deceleration of the motor.

**Parameters** `rpm_minute` – revolution per minute per minute

`set_max_rpm(rpm: int) → None`

Set the maximum RPM.

**Parameters** `rpm` – revolution per minute ( $0 < rpm \leq \text{RPM\_MAX}$ )

**Raises** `ILS2TException` – if RPM is out of range

`set_ramp_type(ramp_type: int = -1) → None`

**Set the ramp type. There are two options available:** 0: linear ramp -1: motor optimized ramp

**Parameters** `ramp_type` – 0: linear ramp | -1: motor optimized ramp

`start() → None`

Start this device.

`stop() → None`

Stop this device. Disables the motor (applies brake), disables access and closes the communication protocol.

`user_steps(steps: int = 16384, revolutions: int = 1) → None`

Define steps per revolution. Default is 16384 steps per revolution. Maximum precision is 32768 steps per revolution.

**Parameters**

- **steps** – number of steps in *revolutions*.
- **revolutions** – number of revolutions corresponding to *steps*.

`write_absolute_position(position: int) → None`

Write instruction to turn the motor until it reaches the absolute position. This function does not enable or disable the motor automatically.

**Parameters** `position` – absolute position of motor in user defined steps.

`write_relative_step(steps: int) → None`

Write instruction to turn the motor the relative amount of steps. This function does not enable or disable the motor automatically.

**Parameters** `steps` – Number of steps to turn the motor.

```
class ILS2TConfig(rpm_max_init: numbers.Integral = 1500, wait_sec_post_enable: Union[int, float] = 1,
                 wait_sec_max_disable: Union[int, float] = 10, wait_sec_post_cannot_disable: Union[int,
                 float] = 1, wait_sec_post_relative_step: Union[int, float] = 2,
                 wait_sec_post_absolute_position: Union[int, float] = 2)
```

Bases: object

Configuration for the ILS2T stepper motor device.

`clean_values()`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**rpm\_max\_init: numbers.Integral = 1500**

initial maximum RPM for the motor, can be set up to 3000 RPM. The user is allowed to set a new max RPM at runtime using *ILS2T.set\_max\_rpm()*, but the value must never exceed this configuration setting.

**wait\_sec\_max\_disable: Union[int, float] = 10**

**wait\_sec\_post\_absolute\_position: Union[int, float] = 2**

**wait\_sec\_post\_cannot\_disable: Union[int, float] = 1**

**wait\_sec\_post\_enable: Union[int, float] = 1**

**wait\_sec\_post\_relative\_step: Union[int, float] = 2**

**exception ILS2TException**

Bases: *hvl\_ccb.dev.base.DeviceException*

Exception to indicate problems with the SE ILS2T stepper motor.

**class ILS2TModbusTcpCommunication(configuration)**

Bases: *hvl\_ccb.comm.modbus\_tcp.ModbusTcpCommunication*

Specific implementation of Modbus/TCP for the Schneider Electric ILS2T stepper motor.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class ILS2TModbusTcpCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], unit: int = 255, port: int = 502)**

Bases: *hvl\_ccb.comm.modbus\_tcp.ModbusTcpCommunicationConfig*

Configuration dataclass for Modbus/TCP communication specific for the Schneider Electric ILS2T stepper motor.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**unit: int = 255**

The unit has to be 255 such that IO scanning mode works.

**class ILS2TRegAddr**(*value*)

Bases: enum.IntEnum

Modbus Register Adresses for for Schneider Electric ILS2T stepper drive.

**ACCESS\_ENABLE = 282**

**FLT\_INFO = 15362**

**FLT\_MEM\_DEL = 15112**

**FLT\_MEM\_RESET = 15114**

**IO\_SCANNING = 6922**

**JOGN\_FAST = 10506**

**JOGN\_SLOW = 10504**

**POSITION = 7706**

**RAMP\_ACC = 1556**

**RAMP\_DECEL = 1558**

**RAMP\_N\_MAX = 1554**

**RAMP\_TYPE = 1574**

**SCALE = 1550**

**TEMP = 7200**

**VOLT = 7198**

**class ILS2TRegDatatype**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: aenum.Enum

Modbus Register Datatypes for Schneider Electric ILS2T stepper drive.

From the manual of the drive:

datatype	byte	min	max
INT8	1 Byte	-128	127
UINT8	1 Byte	0	255
INT16	2 Byte	-32_768	32_767
UINT16	2 Byte	0	65_535
INT32	4 Byte	-2_147_483_648	2_147_483_647
UINT32	4 Byte	0	4_294_967_295
BITS	just 32bits	N/A	N/A

**INT32 = (-2147483648, 2147483647)**

**is\_in\_range**(value: int) → bool

**exception IoScanningModeValueError**

Bases: *hvl\_ccb.dev.se\_ils2t.ILS2TException*


Exception to indicate that the selected IO scanning mode is invalid.

**exception ScalingFactorValueError**

Bases: *hvl\_ccb.dev.se\_ils2t.ILS2TException*

Exception to indicate that a scaling factor value is invalid.

**hvl\_ccb.dev.sst\_luminox**

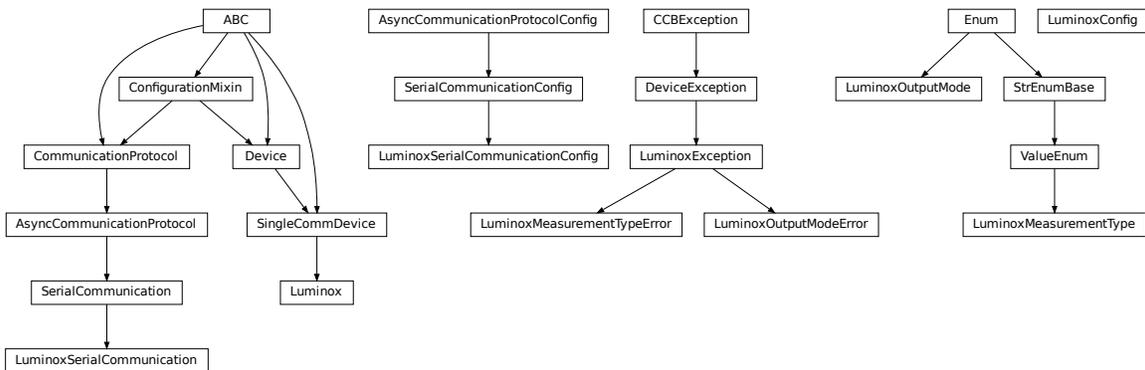

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.

Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

**class Luminox**(com, dev\_config=None)

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Luminox oxygen sensor device class.

**activate\_output**(mode: *hvl\_ccb.dev.sst\_luminox.LuminoxOutputMode*) → None

activate the selected output mode of the Luminox Sensor. :param mode: polling or streaming

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**query\_polling**(*measurement: Union[str, hvl\_ccb.dev.sst\_luminox.LuminoxMeasurementType]*) → Union[Dict[Union[str, *hvl\_ccb.dev.sst\_luminox.LuminoxMeasurementType*], Union[float, int, str]], float, int, str]

Query a value or values of Luminox measurements in the polling mode, according to a given measurement type.

**Parameters** **measurement** – type of measurement

**Returns** value of requested measurement

**Raises**

- **ValueError** – when a wrong key for `LuminoxMeasurementType` is provided
- **LuminoxOutputModeError** – when polling mode is not activated
- **LuminoxMeasurementTypeError** – when expected measurement value is not read

**read\_streaming**() → Dict[Union[str, *hvl\_ccb.dev.sst\_luminox.LuminoxMeasurementType*], Union[float, int, str]]

Read values of Luminox in the streaming mode. Convert the single string into separate values.

**Returns** dictionary with `LuminoxMeasurementType.all_measurements_types()` keys and accordingly type-parsed values.

**Raises**

- **LuminoxOutputModeError** – when streaming mode is not activated
- **LuminoxMeasurementTypeError** – when any of expected measurement values is not read

**start**() → *None*

Start this device. Opens the communication protocol.

**stop**() → *None*

Stop the device. Closes also the communication protocol.

**class LuminoxConfig**(*wait\_sec\_post\_activate: Union[int, float] = 0.5, wait\_sec\_trials\_activate: Union[int, float] = 0.1, nr\_trials\_activate: int = 5*)

Bases: object

Configuration for the SST Luminox oxygen sensor.

**clean\_values**()

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**nr\_trials\_activate: int = 5**

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**wait\_sec\_post\_activate: Union[int, float] = 0.5**

**wait\_sec\_trials\_activate: Union[int, float] = 0.1**

**exception LuminoxException**

Bases: [hvl\\_ccb.dev.base.DeviceException](#)

General Exception for Luminox Device.

**class LuminoxMeasurementType**(*value=<no\_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Measurement types for *LuminoxOutputMode.polling*.

The *all\_measurements* type will read values for the actual measurement types as given in *LuminoxOutputMode.all\_measurements\_types()*; it parses multiple single values using regexp's for other measurement types, therefore, no regexp is defined for this measurement type.

**all\_measurements = 'A'**

**classmethod all\_measurements\_types()** → Tuple[[hvl\\_ccb.dev.sst\\_luminox.LuminoxMeasurementType](#), ...]

A tuple of *LuminoxMeasurementType* enum instances which are actual measurements, i.e. not date of manufacture or software revision.

**barometric\_pressure = 'P'**

**property command: str**

**date\_of\_manufacture = '# 0'**

**parse\_read\_measurement\_value**(*read\_txt: str*) → Union[Dict[Union[str, [hvl\\_ccb.dev.sst\\_luminox.LuminoxMeasurementType](#)], Union[float, int, str]], float, int, str]

**partial\_pressure\_o2 = 'O'**

**percent\_o2 = '%'**

**sensor\_status = 'e'**

**serial\_number = '# 1'**

```
software_revision = '# 2'
```

```
temperature_sensor = 'T'
```

**LuminosMeasurementTypeDict**

A typing hint for a dictionary holding `LuminosMeasurementType` values. Keys are allowed as strings because `LuminosMeasurementType` is of a `StrEnumBase` type.

alias of `Dict[Union[str, LuminosMeasurementType], Union[float, int, str]]`

**exception LuminosMeasurementTypeError**

Bases: `hvl_ccb.dev.sst_luminos.LuminosException`

Wrong measurement type for requested data

**LuminosMeasurementTypeValue**

A typing hint for all possible `LuminosMeasurementType` values as read in either streaming mode or in a polling mode with `LuminosMeasurementType.all_measurements`.

Beware: has to be manually kept in sync with `LuminosMeasurementType` instances `cast_type` attribute values.

alias of `Union[float, int, str]`

**class LuminosOutputMode(value)**

Bases: `enum.Enum`

output mode.

```
polling = 1
```

```
streaming = 0
```

**exception LuminosOutputModeError**

Bases: `hvl_ccb.dev.sst_luminos.LuminosException`

Wrong output mode for requested data

**class LuminosSerialCommunication(configuration)**

Bases: `hvl_ccb.comm.serial.SerialCommunication`

Specific communication protocol implementation for the SST Luminos oxygen sensor. Already predefines device-specific protocol parameters in config.

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class LuminosSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',  
                                         encoding_error_handling: str = 'strict',  
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,  
                                         default_n_attempts_read_text_nonempty: int = 10, port:  
                                         Union[str, NoneType] = None, baudrate: int = 9600, parity:  
                                         Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =  
                                         <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,  
                                         hvl_ccb.comm.serial.SerialCommunicationStopbits] =  
                                         <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,  
                                         hvl_ccb.comm.serial.SerialCommunicationBytesize] =  
                                         <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:  
                                         Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

```
baudrate: int = 9600
```

Baudrate for SST Luminos is 9600 baud

**bytesize:** `Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8`

One byte is eight bits long

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** `Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'`

SST Luminos does not use parity

**classmethod required\_keys**() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** `Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1`

SST Luminos does use one stop bit

**terminator:** `bytes = b'\r\n'`

The terminator is CR LF

**timeout:** `Union[int, float] = 3`

use 3 seconds timeout as default

## [hvl\\_ccb.dev.utils](#)

Poller

**class Poller**(*spoll\_handler*: Callable, *polling\_delay\_sec*: Union[int, float] = 0, *polling\_interval\_sec*: Union[int, float] = 1, *polling\_timeout\_sec*: Optional[Union[int, float]] = None)

Bases: object

Poller class wrapping `concurrent.futures.ThreadPoolExecutor` which enables passing of results and errors out of the polling thread.

**is\_polling()** → bool

Check if device status is being polled.

**Returns** *True* when polling thread is set and alive

**start\_polling()** → bool

Start polling.

**Returns** *True* if was not polling before, *False* otherwise

**stop\_polling()** → bool

Stop polling.

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

**Returns** *True* if was polling before, *False* otherwise, and last result of the polling function call.

**Raises** polling function exceptions


**wait\_for\_polling\_result()**

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

**Returns** polling function result

**Raises** polling function errors

#### hvl\_ccb.dev.visa

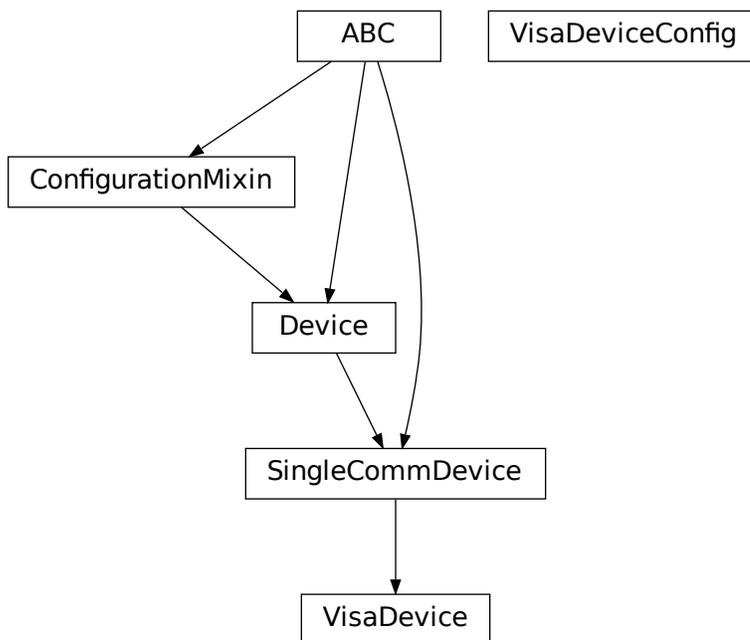

```
class VisaDevice(com: Union[hvl_ccb.comm.visa.VisaCommunication,
                 hvl_ccb.comm.visa.VisaCommunicationConfig, dict], dev_config:
                 Optional[Union[hvl_ccb.dev.visa.VisaDeviceConfig, dict]] = None)
```

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Device communicating over the VISA protocol using VisaCommunication.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()** → Type[*hvl\_ccb.comm.visa.VisaCommunication*]

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns** the VisaCommunication class

**get\_error\_queue()** → str

Read out error queue and logs the error.

**Returns** Error string

**get\_identification()** → str

Queries “\*IDN?” and returns the identification string of the connected device.

**Returns** the identification string of the connected device

**reset()** → *None*

Send “\*RST” and “\*CLS” to the device. Typically sets a defined state.

**spoll\_handler()**

Reads the status byte and decodes it. The status byte STB is defined in IEEE 488.2. It provides a rough overview of the instrument status.

**Returns**

**start()** → *None*

Start the VisaDevice. Sets up the status poller and starts it.

**Returns**

**stop()** → *None*

Stop the VisaDevice. Stops the polling thread and closes the communication protocol.

**Returns**

**wait\_operation\_complete**(*timeout: Optional[float] = None*) → bool

Waits for a operation complete event. Returns after timeout [s] has expired or the operation complete event has been caught.

**Parameters timeout** – Time in seconds to wait for the event; *None* for no timeout.

**Returns** True, if OPC event is caught, False if timeout expired

```
class VisaDeviceConfig(spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2)
```

Bases: *hvl\_ccb.dev.visa.\_VisaDeviceConfigDefaultsBase, hvl\_ccb.dev.visa.*

*\_VisaDeviceConfigBase*

Configdataclass for a VISA device.

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


**Module contents**

Devices subpackage.

**hvl\_ccb.utils**

**Submodules**

**hvl\_ccb.utils.conversion\_sensor**

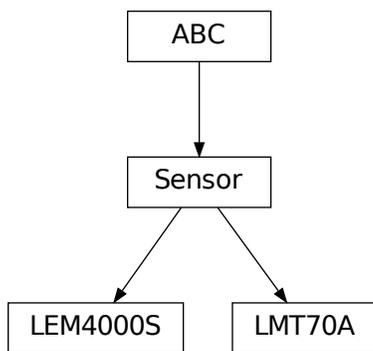

Sensors that are used by the devices implemented in the CCB

**class LEM4000S**(*shunt: float = 1.2, calibration\_factor: float = 1*)  
Bases: [hvl\\_ccb.utils.conversion\\_sensor.Sensor](#)

Converts the output voltage (V) to the measured current (A) when using a LEM Current transducer LT 4000-S

```
CONVERSION: ClassVar[int] = 5000
```

```
calibration_factor: float = 1
```

```
convert(value, **kwargs)
```

```
shunt: float = 1.2
```

```
class LMT70A(temperature_unit: hvl_ccb.utils.conversion_unit.Temperature = Temperature.C)
```

```
Bases: hvl_ccb.utils.conversion_sensor.Sensor
```

Converts the output voltage (V) to the measured temperature (default °C) when using a TI Precision Analog Temperature Sensor LMT70(A)

```
LUT: ClassVar[numpy.ndarray[Any,
numpy.dtype[numpy.typing._generic_alias.ScalarType]]] = array([[ -55. , 1.375219],
[ -50. , 1.350441], [ -40. , 1.300593], [ -30. , 1.250398], [ -20. , 1.199884], [ -10. ,
1.14907 ], [  0. , 1.097987], [ 10. , 1.046647], [ 20. , 0.99505 ], [ 30. ,
0.943227], [ 40. , 0.891178], [ 50. , 0.838882], [ 60. , 0.78636 ], [ 70. ,
0.733608], [ 80. , 0.680654], [ 90. , 0.62749 ], [100. , 0.574117], [110. ,
0.520551], [120. , 0.46676 ], [130. , 0.412739], [140. , 0.358164], [150. ,
0.302785]])
```

```
convert(value, **kwargs)
```

```
temperature_unit: hvl_ccb.utils.conversion_unit.Temperature = 'C'
```


```
class Sensor
```

```
Bases: abc.ABC
```

The BaseClass ‘Sensor’ is designed as a parent for all Sensors. Each attribute must be added to ‘\_\_setattr\_\_’, so that the value is verified each time the value is changed. It is important to mark attributes that should be constant with ‘typing.ClassVar[...]’. Together with ‘super().\_\_setattr\_\_(name, value)’, this guarantees that the values are protected and cannot be altered by the user.

```
abstract convert(value, **kwargs)
```

### hvl\_ccb.utils.conversion\_unit

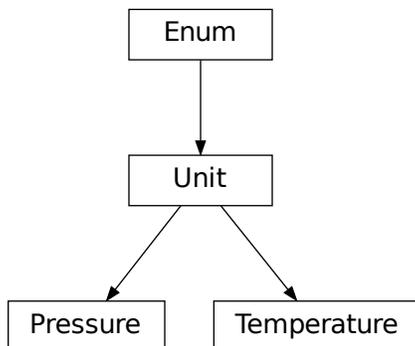

example Kelvin <-> Celsius

Unit conversion, within in the same group of units, for

**class Pressure**(*value*)

Bases: *hvl\_ccb.utils.conversion\_unit.Unit*

An enumeration.

**ATM** = 'atm'

**ATMOSPHERE** = 'atm'

**BAR** = 'bar'

**MILLIMETER\_MERCURY** = 'mmHg'

**MMHG** = 'mmHg'

**PA** = 'Pa'

**PASCAL** = 'Pa'

**POUNDS\_PER\_SQUARE\_INCH** = 'psi'

**PSI** = 'psi'

**TORR** = 'torr'

**class Temperature**(*value*)

Bases: *hvl\_ccb.utils.conversion\_unit.Unit*

An enumeration.

**C** = 'C'

**CELSIUS** = 'C'

**F** = 'F'

**FAHRENHEIT** = 'F'

**K** = 'K'

**KELVIN** = 'K'

**class Unit**(*value*)

Bases: `enum.Enum`


An enumeration.

**abstract classmethod convert**(*value*, *\*\*kwargs*)

**preserve\_type**(*func*)

This wrapper preserves the first order type of the input. Upto now the type of the data stored in a list, tuple, array or dict is not preserved. Integer will be converted to float!

## hvl\_ccb.utils.enum

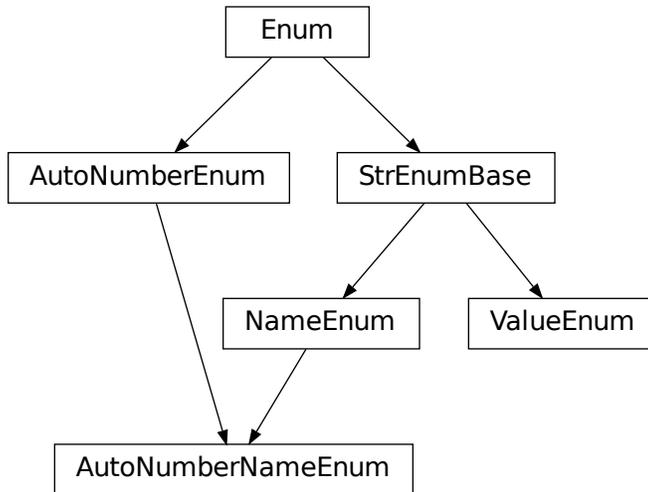

**class AutoNumberNameEnum**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.NameEnum`, `aenum.AutoNumberEnum`

Auto-numbered enum with names used as string representation, and with lookup and equality based on this representation.

**class NameEnum**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.StrEnumBase`

Enum with names used as string representation, and with lookup and equality based on this representation.

**class StrEnumBase**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `aenum.Enum`

String representation-based equality and lookup.

**class ValueEnum**(*value=<no\_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.StrEnumBase`

Enum with string representation of values used as string representation, and with lookup and equality based on this representation.

Attention: to avoid errors, best use together with *unique* enum decorator.

## hvl\_ccb.utils.typing

Additional Python typing module utilities

### ConvertibleTypes

Typing hint for data type that can be used in conversion

alias of `Union[int, float, List[Union[int, float]], Tuple[Union[int, float]], Dict[str, Union[int, float]], numpy.ndarray]`

### Number

Typing hint auxiliary for a Python base number types: *int* or *float*.

alias of `Union[int, float]`

### check\_generic\_type(value, type\_, name='instance')

Check if *value* is of a generic type *type\_*. Raises *TypeError* if it's not.

#### Parameters

- **name** – name to report in case of an error
- **value** – value to check
- **type** – generic type to check against

### is\_generic\_type\_hint(type\_)

Check if class is a generic type, for example *Union[int, float]*, *List[int]*, or *List*.

**Parameters** **type** – type to check

## hvl\_ccb.utils.validation

**validate\_and\_resolve\_host**(*host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]*, *logger: Optional[logging.Logger] = None*) → *str*

**validate\_bool**(*x\_name: str*, *x: object*, *logger: Optional[logging.Logger] = None*) → *None*  
Validate if given input *x* is a *bool*.

#### Parameters

- **x\_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as boolean
- **logger** – logger of the calling submodule

**Raises** **TypeError** – when the validated input does not have boolean type

**validate\_number**(*x\_name: str*, *x: object*, *limits: typing.Optional[typing.Tuple] = (None, None)*, *number\_type: typing.Union[typing.Type[typing.Union[int, float]], typing.Tuple[typing.Type[typing.Union[int, float]], ...]] = (<class 'int'>, <class 'float'>)*, *logger: typing.Optional[logging.Logger] = None*) → *None*

Validate if given input *x* is a number of given *number\_type* type, with value between given *limits[0]* and *limits[1]* (inclusive), if not *None*. For array-like objects (*npt.NDArray*, *List*, *Tuple*, *Dict*) it is checked if all elements are within the limits and have the correct type.

#### Parameters

- **x\_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as number of given type within given range
- **logger** – logger of the calling submodule

- **limits** – [lower, upper] limit, with *None* denoting no limit: [-inf, +inf]
- **number\_type** – expected type or tuple of types of a number, by default (*int, float*)

#### Raises


- **TypeError** – when the validated input does not have expected type
- **ValueError** – when the validated input has correct number type but is not within given range

`validate_tcp_port(port: int, logger: Optional[logging.Logger] = None)`

## Module contents

### 4.1.2 Submodules

#### hvl\_ccb.configuration

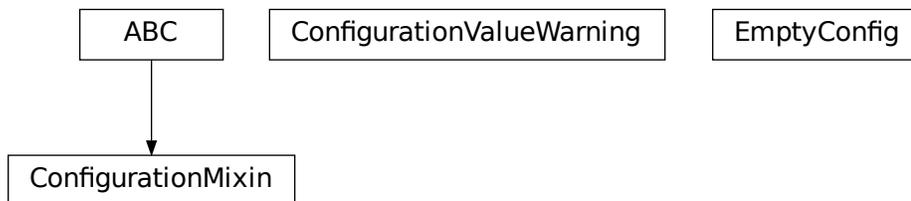

#### Facilities

providing classes for handling configuration for communication protocols and devices.

**class ConfigurationMixin**(*configuration*)

Bases: abc.ABC

Mixin providing configuration to a class.

**property config**

ConfigDataclass property.

**Returns** the configuration

**abstract static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**configuration\_save\_json**(*path: str*) → *None*

Save current configuration as JSON file.

**Parameters path** – path to the JSON file.

**classmethod from\_json**(*filename: str*)

Instantiate communication protocol using configuration from a JSON file.

**Parameters filename** – Path and filename to the JSON configuration

**exception ConfigurationValueWarning**

Bases: UserWarning

User warnings category for values of *@configdataclass* fields.**class EmptyConfig**

Bases: object

Empty configuration dataclass.

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True****classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


**Returns** a list of strings containing all required keys.**configdataclass(*direct\_decoration=None*, *frozen=True*)**Decorator to make a class a configdataclass. Types in these dataclasses are enforced. Implement a function *clean\_values(self)* to do additional checking on value ranges etc.It is possible to inherit from a configdataclass and re-decorate it with *@configdataclass*. In a subclass, default values can be added to existing fields. Note: adding additional non-default fields is prone to errors, since the order has to be respected through the whole chain (first non-default fields, only then default-fields).**Parameters** **frozen** – defaults to True. False allows to later change configuration values. Attention: if configdataclass is not frozen and a value is changed, typing is not enforced anymore!

**hvl\_ccb.exception**

CCBException

Introduce a common base exception for the CCB

**exception CCBException**  
Bases: Exception

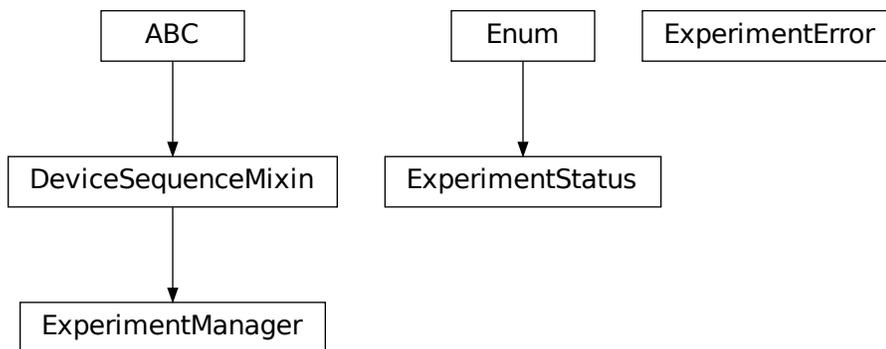**hvl\_ccb.experiment\_manager**

Main module containing the top level ExperimentManager class. Inherit from this class to implement your own experiment functionality in another project and it will help you start, stop and manage your devices.

**exception ExperimentError**  
Bases: Exception

Exception to indicate that the current status of the experiment manager is on ERROR and thus no operations can be made until reset.

**class ExperimentManager(\*args, \*\*kwargs)**  
Bases: *hvl\_ccb.dev.base.DeviceSequenceMixin*

Experiment Manager can start and stop communication protocols and devices. It provides methods to queue commands to devices and collect results.

**add\_device(name: str, device: hvl\_ccb.dev.base.Device) → None**

Add a new device to the manager. If the experiment is running, automatically start the device. If a device with this name already exists, raise an exception.

**Parameters**

- **name** – is the name of the device.

- **device** – is the instantiated Device object.

Raises *DeviceExistingException* –

**devices\_failed\_start:** Dict[str, *hvl\_ccb.dev.base.Device*]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices\_failed\_stop*.

**devices\_failed\_stop:** Dict[str, *hvl\_ccb.dev.base.Device*]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices\_failed\_start*.

**finish()** → *None*

Stop experimental setup, stop all devices.

**is\_error()** → bool

Returns true, if the status of the experiment manager is *error*.

**Returns** True if on error, false otherwise

**is\_finished()** → bool

Returns true, if the status of the experiment manager is *finished*.

**Returns** True if finished, false otherwise

**is\_running()** → bool

Returns true, if the status of the experiment manager is *running*.

**Returns** True if running, false otherwise

**run()** → *None*

Start experimental setup, start all devices.

**start()** → *None*

Alias for ExperimentManager.run()

**property status:** *hvl\_ccb.experiment\_manager.ExperimentStatus*

Get experiment status.

**Returns** experiment status enum code.

**stop()** → *None*

Alias for ExperimentManager.finish()

**class ExperimentStatus**(*value*)

Bases: enum.Enum

Enumeration for the experiment status

**ERROR** = 5

**FINISHED** = 4

**FINISHING** = 3

**INITIALIZED** = 0

**INITIALIZING** = -1

**RUNNING** = 2

**STARTING** = 1

### 4.1.3 Module contents

Top-level package for HVL Common Code Base.



## CONTRIBUTING

Contributions are welcome, and they are greatly appreciated! Every little bit helps, and credit will always be given.

You can contribute in many ways:

### 5.1 Types of Contributions

#### 5.1.1 Report Bugs

Report bugs at [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/issues](https://gitlab.com/ethz_hvl/hvl_ccb/issues).

If you are reporting a bug, please include:

- Your operating system name and version.
- Any details about your local setup that might be helpful in troubleshooting.
- Detailed steps to reproduce the bug.

#### 5.1.2 Fix Bugs

Look through the GitLab issues for bugs. Anything tagged with “bug” and “help wanted” is open to whoever wants to implement it.

#### 5.1.3 Implement Features

Look through the GitLab issues for features. Anything tagged with “enhancement” and “help wanted” is open to whoever wants to implement it.

#### 5.1.4 Write Documentation

HVL Common Code Base could always use more documentation, whether as part of the official HVL Common Code Base docs, in docstrings, or even on the web in blog posts, articles, and such.

## 5.1.5 Submit Feedback

The best way to send feedback is to file an issue at [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/issues](https://gitlab.com/ethz_hvl/hvl_ccb/issues).

If you are proposing a feature:

- Explain in detail how it would work.
- Keep the scope as narrow as possible, to make it easier to implement.
- Remember that this is a volunteer-driven project, and that contributions are welcome :)

## 5.2 Get Started!

Ready to contribute? Here's how to set up *hvl\_ccb* for local development.

1. Clone *hvl\_ccb* repo from GitLab.

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

2. Install your local copy into a virtualenv. Assuming you have *virtualenvwrapper* installed, this is how you set up your fork for local development:

```
$ mkvirtualenv hvl_ccb
$ cd hvl_ccb/
$ pip install -e .[all]
$ pip install -r requirements_dev.txt
```

3. Create a branch for local development:

```
$ git checkout -b name-of-your-bugfix-or-feature
```

Now you can make your changes locally.

4. When you're done making changes, check that your changes pass *flake8* and the tests, including testing other Python versions with *tox*:

```
$ flake8 hvl_ccb tests
$ python setup.py test or py.test
$ tox
```

To get *flake8* and *tox*, just *pip* install them into your virtualenv. You can also use the provided *make*-like shell script to run *flake8* and tests:

```
$ ./make.sh lint
$ ./make.sh test
```

5. Commit your changes and push your branch to GitLab:

```
$ git add .
$ git commit -m "Your detailed description of your changes."
$ git push origin name-of-your-bugfix-or-feature
```

6. Submit a merge request through the GitLab website.

## 5.3 Merge Request Guidelines

Before you submit a merge request, check that it meets these guidelines:

1. The merge request should include tests.
2. If the merge request adds functionality, the docs should be updated. Put your new functionality into a function with a docstring, and add the feature to the list in README.rst.
3. The merge request should work for Python 3.7. Check [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/merge\\_requests](https://gitlab.com/ethz_hvl/hvl_ccb/merge_requests) and make sure that the tests pass for all supported Python versions.

## 5.4 Tips

- To run tests from a single file:

```
$ py.test tests/test_hvl_ccb.py
```

or a single test function:

```
$ py.test tests/test_hvl_ccb.py::test_command_line_interface
```

- If your tests are slow, profile them using the pytest-profiling plugin:

```
$ py.test tests/test_hvl_ccb.py --profile
```

or for a graphical overview (you need a SVG image viewer):

```
$ py.test tests/test_hvl_ccb.py --profile-svg
$ open prof/combined.svg
```

- To add dependency, edit appropriate `*requirements` variable in the `setup.py` file and re-run:

```
$ python setup.py develop
```

- To generate a PDF version of the Sphinx documentation instead of HTML use:

```
$ rm -rf docs/hvl_ccb.rst docs/modules.rst docs/_build && sphinx-apidoc -o docs/hvl_
↪ccb && python -msphinx -M latexpdf docs/ docs/_build
```

This command can also be run through the make-like shell script:

```
$ ./make.sh docs-pdf
```

This requires a local installation of a LaTeX distribution, e.g. MikTeX.

## 5.5 Deploying

A reminder for the maintainers on how to deploy. Create `release-N.M.K` branch. Make sure all your changes are committed. Update or create entry in `HISTORY.rst` file, and, if applicable, update `AUTHORS.rst` file, update features tables in `README.rst` file, and update API docs:

```
$ make docs
```

Commit all of the above, except for the `docs/hvl_ccb.dev.picotech_pt104.rst`, and then run:

```
$ bumpversion patch # possible: major / minor / patch
$ git push
$ git push --tags
```

Go to <https://readthedocs.org/projects/hvl-ccb/builds/> and check if RTD docs build for the pushed tag passed.

Wait for the CI pipeline to finish successfully; afterwards, run a release check:

```
$ make release-check
```

Finally, prepare and push a release:

```
$ make release
```

Merge the release branch into master and devel branches with `--no-ff` flag and delete the release branch:

```
$ git switch master
$ git merge --no-ff release-N.M.K
$ git push
$ git switch devel
$ git merge --no-ff release-N.M.K
$ git push
$ git push --delete origin release-N.M.K
$ git branch --delete release-N.M.K
```

Finally, prepare GitLab release and cleanup the corresponding milestone:

1. go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/-/tags/](https://gitlab.com/ethz_hvl/hvl_ccb/-/tags/), select the latest release tag, press “Edit release notes” and add the release notes (copy a corresponding entry from `HISTORY.rst` file with formatting adjusted from ReStructuredText to Markdown); press “Save changes”;
2. go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/-/releases](https://gitlab.com/ethz_hvl/hvl_ccb/-/releases), select the latest release, press “Edit this release” and under “Milestones” select the corresponding milestone; press “Save changes”;
3. go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/-/milestones](https://gitlab.com/ethz_hvl/hvl_ccb/-/milestones), make sure that it is 100% complete (otherwise, create a next patch-level milestone and assign it to the ongoing Issues and Merge Requests therein); press “Close Milestone”.

## 6.1 Maintainers

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>

## 6.2 Authors

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- David Graber <dev@davidgraber.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Alise Chachereau <chachereau@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- David Taylor <dtaylor@ethz.ch>
- Joseph Engelbrecht <engelbrecht@eeh.ee.ethz.ch>

## 6.3 Contributors

- Luca Nembrini <lucane@student.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>
- Raphael Faerber <raphael.ferber@eeh.ee.ethz.ch>
- Ruben Stadler <rstadler@student.ethz.ch>
- Hanut Vemulapalli <vemulapalli@eeh.ee.ethz.ch>



## HISTORY

### 7.1 0.10.0 (2022-01-17)

- Reimplementation of the Cube (before known as Supercube)
- **new names:**
  - Supercube Typ B -> BaseCube
  - Supercube Typ A -> PICube (power inverter Cube)
- **new import:**
  - `from hvl_ccb.dev.supercube import SupercubeB -> from hvl_ccb.dev.cube import BaseCube`
- **new programming style:**
  - getter / setter methods -> properties
  - e.g. get: `cube.get_support_output(port=1, contact=1) -> cube.support_1.output_1`
  - e.g. set: `cube.get_support_output(port=1, contact=1, state=True) -> cube.support_1.output_1 = True`
- unify Exceptions of Cube
- implement Fast Switch-Off of Cube
- remove method `support_output_impulse`
- all active alarms can now be queried `cube.active_alarms()`
- alarms will now result in different logging levels depending on the seriousness of the alarm.
- introduction of limits for slope and safety limit for RedReady
- during the startup the CCB will update the time of the cube.
- verification of inputs
- polarity of DC voltage
- Switch from `python-opcua` to `opcua-asyncio` (former package is no longer maintained)

## 7.2 0.9.0 (2022-01-07)

- New device: Highland T560 digital delay and pulse generator over Telnet.
- **Rework of the Technix Capacitor Charger.**
  - Moved into a separate sub-package
  - NEW import over `import hvl_ccb.dev.technix as XXX`
  - Slightly adapted behaviour
- Add `validate_tcp_port` to validate port number.
- **Add `validate_and_resolve_host` to validate and resolve host names and IPs.**
  - Remove requirement `IPy`
- Add a unified CCB Exception schema for all devices and communication protocols.
- Add data conversion functions to README.
- Update CI and devel images from Debian 10 buster to Debian 11 bullseye.
- Fix typing due to numpy update.
- Fix incorrect overloading of `clean_values()` in classes of type `XCommunicationConfig`.

## 7.3 0.8.5 (2021-11-05)

- Added arbitrary waveform for TiePie signal generation, configurable via `dev.tiepie.generator.TiePieGeneratorConfig.waveform` property.
- In `utils.conversion_sensor`: improvements for class constants; removed SciPy dependency.
- Added Python 3.10 support.

## 7.4 0.8.4 (2021-10-22)

- `utils.validation.validate_number` extension to handle NumPy arrays and array-like objects.
- `utils.conversion_unit` utility classes handle correctly `NamedTuple` instances.
- `utils.conversion_sensor` and `utils.conversion_unit` code simplification (no `transfer_function_order` attribute) and cleanups.
- Fixed incorrect error logging in `configuration.configdataclass`.
- `comm.telnet.TelnetCommunication` tests fixes for local run errors.

## 7.5 0.8.3 (2021-09-27)

- New data conversion functions in `utils.conversion_sensor` and `utils.conversion_unit` modules. Note: to use these functions you must install `hvl_ccb` with extra requirement, either `hvl_ccb[conversion]` or `hvl_ccb[all]`.
- Improved documentation with respect to installation of external libraries.

## 7.6 0.8.2 (2021-08-27)

- **New functionality in `dev.labjack.LabJack`:**
  - configure clock and send timed pulse sequences
  - set DAC/analog output voltage
- Bugfix: ignore random bits sent by to `dev.newport.NewportSMC100PP` controller during start-up/powering-up.

## 7.7 0.8.1 (2021-08-13)

- Add Python version check (min version error; max version warning).
- Daily checks for upstream dependencies compatibility and devel environment improvements.

## 7.8 0.8.0 (2021-07-02)

- TCP communication protocol.
- Lauda PRO RP 245 E circulation thermostat device over TCP.
- Pico Technology PT-104 Platinum Resistance Data Logger device as a wrapper of the Python bindings for the PicoSDK.
- In `com.visa.VisaCommunication`: periodic status polling when VISA/TCP keep alive connection is not supported by a host.

## 7.9 0.7.1 (2021-06-04)

- New `utils.validation` submodule with `validate_bool` and `validate_number` utilities extracted from internal use within a `dev.tiepie` subpackage.
- **In `comm.serial.SerialCommunication`:**
  - strict encoding errors handling strategy for subclasses,
  - user warning for a low communication timeout value.

## 7.10 0.7.0 (2021-05-25)

- The `dev.tiepie` module was splitted into a subpackage with, in particular, submodules for each of the device types – `oscilloscope`, `generator`, and `i2c` – and with backward-incompatible direct imports from the submodules.
- **In `dev.technix`:**
  - fixed communication crash on nested status byte query;
  - added enums for GET and SET register commands.
- Further minor logging improvements: added missing module level logger and removed some error logs in `except` blocks used for a flow control.
- In `examples/` folder renamed consistently all the examples.
- In API documentation: fix incorrect links mapping on inheritance diagrams.

## 7.11 0.6.1 (2021-05-08)

- **In `dev.tiepie`:**
  - dynamically set `oscilloscope`'s channel limits in `OscilloscopeChannelParameterLimits`: `input_range` and `trigger_level_abs`, incl. update of latter on each change of `input_range` value of a `TiePieOscilloscopeChannelConfig` instances;
  - quick fix for opening of combined instruments by disabling `OscilloscopeParameterLimits`. `trigger_delay` (an advanced feature);
  - enable automatic devices detection to be able to find network devices with `TiePieOscilloscope`. `list_devices()`.
- Fix `examples/example_labjack.py`.
- Improved logging: consistently use module level loggers, and always log exception tracebacks.
- Improve API documentation: separate pages per modules, each with an inheritance diagram as an overview.

## 7.12 0.6.0 (2021-04-23)

- Technix capacitor charger using either serial connection or Telnet protocol.
- **Extensions, improvements and fixes in existing devices:**
  - **In `dev.tiepie.TiePieOscilloscope`:**
    - \* redesigned measurement start and data collection API, incl. time out argument, with no/infinite time out option;
    - \* trigger allows now a no/infinite time out;
    - \* record length and trigger level were fixed to accept, respectively, floating point and integer numbers;
    - \* fixed resolution validation bug;
  - `dev.heinzinger.HeinzingerDI` and `dev.rs_rto1024.RTO1024` instances are now resilient to multiple `stop()` calls.

- In `dev.crylas.CryLasLaser`: default configuration timeout and polling period were adjusted;
- Fixed PSI9080 example script.
- **Package and source code improvements:**
  - Update to backward-incompatible `pyvisa-py>=0.5.2`. Developers, do update your local development environments!
  - External libraries, like LibTiePie SDK or LJM Library, are now not installed by default; they are now extra installation options.
  - Added Python 3.9 support.
  - Improved number formatting in logs.
  - Typing improvements and fixes for `mypy>=0.800`.

## 7.13 0.5.0 (2020-11-11)

- TiePie USB oscilloscope, generator and I2C host devices, as a wrapper of the Python bindings for the LibTiePie SDK.
- a FuG Elektronik Power Supply (e.g. Capacitor Charger HCK) using the built-in ADDAT controller with the Probus V protocol over a serial connection
- All devices polling status or measurements use now a `dev.utils.Poller` utility class.
- **Extensions and improvements in existing devices:**
  - In `dev.rs_rto1024.RT01024`: added Channel state, scale, range, position and offset accessors, and measurements activation and read methods.
  - In `dev.sst_luminox.Luminox`: added querying for all measurements in polling mode, and made output mode activation more robust.
  - In `dev.newport.NewportSMC100PP`: an error-prone `wait_until_move_finished` method of replaced by a fixed waiting time, device operations are now robust to a power supply cut, and device restart is not required to apply a start configuration.
- **Other minor improvements:**
  - Single failure-safe starting and stopping of devices sequenced via `dev.base.DeviceSequenceMixin`.
  - Moved `read_text_nonempty` up to `comm.serial.SerialCommunication`.
  - Added development Dockerfile.
  - Updated package and development dependencies: `pymodbus`, `pytest-mock`.

## 7.14 0.4.0 (2020-07-16)

- **Significantly improved new Supercube device controller:**
  - more robust error-handling,
  - status polling with generic Poller helper,
  - messages and status boards.
  - tested with a physical device,
- Improved OPC UA client wrapper, with better error handling, incl. re-tries on `concurrent.futures.TimeoutError`.
- SST Luminox Oxygen sensor device controller.
- **Backward-incompatible changes:**
  - `CommunicationProtocol.access_lock` has changed type from `threading.Lock` to `threading.RLock`.
  - `ILS2T.relative_step` and `ILS2T.absolute_position` are now called, respectively, `ILS2T.write_relative_step` and `ILS2T.write_absolute_position`.
- **Minor bugfixes and improvements:**
  - fix use of max resolution in `Labjack.set_ain_resolution()`,
  - resolve ILS2T devices relative and absolute position setters race condition,
  - added acoustic horn function in the 2015 Supercube.
- **Toolchain changes:**
  - add Python 3.8 support,
  - drop pytest-runner support,
  - ensure compatibility with `labjack_ljm` 2019 version library.

## 7.15 0.3.5 (2020-02-18)

- Fix issue with reading integers from LabJack LJM Library (device's product ID, serial number etc.)
- Fix development requirements specification (tox version).

## 7.16 0.3.4 (2019-12-20)

- **New devices using serial connection:**
  - Heinzinger Digital Interface I/II and a Heinzinger PNC power supply
  - Q-switched Pulsed Laser and a laser attenuator from CryLas
  - Newport SMC100PP single axis motion controller for 2-phase stepper motors
  - Pfeiffer TPG controller (TPG 25x, TPG 26x and TPG 36x) for Compact pressure Gauges
- PEP 561 compatibility and related corrections for static type checking (now in CI)
- **Refactorings:**

- Protected non-thread safe read and write in communication protocols
- Device sequence mixin: start/stop, add/rm and lookup
- *.format()* to f-strings
- more enumerations and a quite some improvements of existing code
- Improved error docstrings (:raises: annotations) and extended tests for errors.

## 7.17 0.3.3 (2019-05-08)

- Use PyPI labjack-ljm (no external dependencies)

## 7.18 0.3.2 (2019-05-08)

- INSTALLATION.rst with LJMPython prerequisite info

## 7.19 0.3.1 (2019-05-02)

- readthedocs.org support

## 7.20 0.3 (2019-05-02)

- Prevent an automatic close of VISA connection when not used.
- Rhode & Schwarz RTO 1024 oscilloscope using VISA interface over `TCP::INSTR`.
- Extended tests incl. messages sent to devices.
- Added Supercube device using an OPC UA client
- Added Supercube 2015 device using an OPC UA client (for interfacing with old system version)

## 7.21 0.2.1 (2019-04-01)

- Fix issue with LJMPython not being installed automatically with setuptools.

## 7.22 0.2.0 (2019-03-31)

- LabJack LJM Library communication wrapper and LabJack device.
- Modbus TCP communication protocol.
- Schneider Electric ILS2T stepper motor drive device.
- Elektro-Automatik PSI9000 current source device and VISA communication wrapper.
- Separate configuration classes for communication protocols and devices.
- Simple experiment manager class.

## 7.23 0.1.0 (2019-02-06)

- Communication protocol base and serial communication implementation.
- Device base and MBW973 implementation.

## INDICES AND TABLES

- `genindex`
- `modindex`
- `search`



## PYTHON MODULE INDEX

### h

hvl\_ccb, 191  
hvl\_ccb.comm, 33  
hvl\_ccb.comm.base, 11  
hvl\_ccb.comm.labjack\_ljm, 15  
hvl\_ccb.comm.modbus\_tcp, 18  
hvl\_ccb.comm.opc, 20  
hvl\_ccb.comm.serial, 23  
hvl\_ccb.comm.tcp, 26  
hvl\_ccb.comm.telnet, 28  
hvl\_ccb.comm.visa, 30  
hvl\_ccb.configuration, 187  
hvl\_ccb.dev, 182  
hvl\_ccb.dev.base, 92  
hvl\_ccb.dev.crylas, 94  
hvl\_ccb.dev.cube, 53  
hvl\_ccb.dev.cube.base, 33  
hvl\_ccb.dev.cube.constants, 38  
hvl\_ccb.dev.cube.picube, 50  
hvl\_ccb.dev.ea\_psi9000, 104  
hvl\_ccb.dev.fug, 108  
hvl\_ccb.dev.heinzinger, 119  
hvl\_ccb.dev.highland\_t560, 58  
hvl\_ccb.dev.highland\_t560.base, 53  
hvl\_ccb.dev.highland\_t560.channel, 55  
hvl\_ccb.dev.highland\_t560.device, 56  
hvl\_ccb.dev.labjack, 125  
hvl\_ccb.dev.lauda, 132  
hvl\_ccb.dev.mbw973, 138  
hvl\_ccb.dev.newport, 142  
hvl\_ccb.dev.pfeiffer\_tpg, 155  
hvl\_ccb.dev.rs\_rto1024, 161  
hvl\_ccb.dev.se\_ils2t, 169  
hvl\_ccb.dev.sst\_luminox, 175  
hvl\_ccb.dev.supercube2015, 73  
hvl\_ccb.dev.supercube2015.base, 58  
hvl\_ccb.dev.supercube2015.constants, 64  
hvl\_ccb.dev.supercube2015.typ\_a, 71  
hvl\_ccb.dev.technix, 79  
hvl\_ccb.dev.technix.base, 74  
hvl\_ccb.dev.technix.device, 77  
hvl\_ccb.dev.tiepie, 91  
hvl\_ccb.dev.tiepie.base, 80  
hvl\_ccb.dev.tiepie.channel, 82  
hvl\_ccb.dev.tiepie.device, 84  
hvl\_ccb.dev.tiepie.generator, 85  
hvl\_ccb.dev.tiepie.i2c, 87  
hvl\_ccb.dev.tiepie.oscilloscope, 88  
hvl\_ccb.dev.tiepie.utils, 90  
hvl\_ccb.dev.utils, 179  
hvl\_ccb.dev.visa, 180  
hvl\_ccb.exception, 189  
hvl\_ccb.experiment\_manager, 189  
hvl\_ccb.utils, 187  
hvl\_ccb.utils.conversion\_sensor, 182  
hvl\_ccb.utils.conversion\_unit, 183  
hvl\_ccb.utils.enum, 185  
hvl\_ccb.utils.typing, 186  
hvl\_ccb.utils.validation, 186



## A

- A (*HeinzingerPNC.UnitCurrent* attribute), 122
- A (*SupercubeOpcEndpoint* attribute), 71
- ABSOLUTE (*TiePieOscilloscopeTriggerLevelMode* attribute), 84
- ABSOLUTE\_POSITION (*ILS2T.ActionsPtp* attribute), 169
- AC (*NewportConfigCommands* attribute), 142
- AC\_100KV (*PowerSetup* attribute), 49
- AC\_150KV (*PowerSetup* attribute), 49
- AC\_200KV (*PowerSetup* attribute), 49
- AC\_50KV (*PowerSetup* attribute), 49
- AC\_DoubleStage\_150kV (*PowerSetup* attribute), 69
- AC\_DoubleStage\_200kV (*PowerSetup* attribute), 69
- AC\_SingleStage\_100kV (*PowerSetup* attribute), 69
- AC\_SingleStage\_50kV (*PowerSetup* attribute), 69
- ACA (*TiePieOscilloscopeChannelCoupling* attribute), 83
- acceleration (*NewportSMC100PPConfig* attribute), 150
- ACCESS\_ENABLE (*ILS2TRegAddr* attribute), 174
- access\_lock (*CommunicationProtocol* attribute), 14
- ACTION\_JOG\_VALUE (*ILS2T* attribute), 169
- activate\_clock\_output() (*T560* method), 56
- activate\_measurements() (*RTO1024* method), 162
- activate\_output() (*Luminos* method), 175
- activated (*BreakdownDetection* attribute), 65
- ACTIVE (*CryLasLaser.AnswersStatus* attribute), 97
- active\_alarms() (*BaseCube* method), 33
- ACTIVE\_HIGH (*Polarity* attribute), 53
- ACTIVE\_LOW (*Polarity* attribute), 54
- actualsetvalue (*FuGProbusVSetRegisters* property), 116
- ACV (*TiePieOscilloscopeChannelCoupling* attribute), 83
- adc\_mode (*FuGProbusVMonitorRegisters* property), 116
- add\_device() (*DeviceSequenceMixin* method), 92
- add\_device() (*ExperimentManager* method), 189
- ADDR\_INCORRECT (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 151
- address (*NewportSMC100PPConfig* attribute), 150
- address (*VisaCommunicationConfig* property), 31
- address() (*VisaCommunicationConfig.InterfaceType* method), 31
- ADMODE (*FuGProbusIVCommands* attribute), 113
- Alarm0 (*AlarmText* attribute), 64
- Alarm1 (*AlarmText* attribute), 64
- Alarm10 (*AlarmText* attribute), 64
- Alarm11 (*AlarmText* attribute), 64
- Alarm12 (*AlarmText* attribute), 64
- Alarm13 (*AlarmText* attribute), 64
- Alarm14 (*AlarmText* attribute), 64
- Alarm17 (*AlarmText* attribute), 64
- Alarm19 (*AlarmText* attribute), 64
- Alarm2 (*AlarmText* attribute), 64
- Alarm20 (*AlarmText* attribute), 64
- Alarm21 (*AlarmText* attribute), 64
- Alarm22 (*AlarmText* attribute), 64
- Alarm3 (*AlarmText* attribute), 64
- Alarm4 (*AlarmText* attribute), 64
- Alarm5 (*AlarmText* attribute), 64
- Alarm6 (*AlarmText* attribute), 64
- Alarm7 (*AlarmText* attribute), 64
- Alarm8 (*AlarmText* attribute), 64
- Alarm9 (*AlarmText* attribute), 64
- ALARM\_1 (*Alarms* attribute), 38
- alarm\_1 (*AlarmsOverview* attribute), 42
- ALARM\_10 (*Alarms* attribute), 38
- alarm\_10 (*AlarmsOverview* attribute), 42
- ALARM\_100 (*Alarms* attribute), 38
- alarm\_100 (*AlarmsOverview* attribute), 42
- ALARM\_101 (*Alarms* attribute), 38
- alarm\_101 (*AlarmsOverview* attribute), 42
- ALARM\_102 (*Alarms* attribute), 38
- alarm\_102 (*AlarmsOverview* attribute), 42
- ALARM\_103 (*Alarms* attribute), 38
- alarm\_103 (*AlarmsOverview* attribute), 42
- ALARM\_104 (*Alarms* attribute), 38
- alarm\_104 (*AlarmsOverview* attribute), 42
- ALARM\_105 (*Alarms* attribute), 38
- alarm\_105 (*AlarmsOverview* attribute), 42
- ALARM\_106 (*Alarms* attribute), 38
- alarm\_106 (*AlarmsOverview* attribute), 43
- ALARM\_107 (*Alarms* attribute), 38
- alarm\_107 (*AlarmsOverview* attribute), 43
- ALARM\_108 (*Alarms* attribute), 38





- alarm\_7 (*AlarmsOverview* attribute), 46
- ALARM\_70 (*Alarms* attribute), 41
- alarm\_70 (*AlarmsOverview* attribute), 46
- ALARM\_71 (*Alarms* attribute), 41
- alarm\_71 (*AlarmsOverview* attribute), 46
- ALARM\_72 (*Alarms* attribute), 41
- alarm\_72 (*AlarmsOverview* attribute), 46
- ALARM\_73 (*Alarms* attribute), 41
- alarm\_73 (*AlarmsOverview* attribute), 46
- ALARM\_74 (*Alarms* attribute), 41
- alarm\_74 (*AlarmsOverview* attribute), 46
- ALARM\_75 (*Alarms* attribute), 41
- alarm\_75 (*AlarmsOverview* attribute), 46
- ALARM\_76 (*Alarms* attribute), 41
- alarm\_76 (*AlarmsOverview* attribute), 46
- ALARM\_77 (*Alarms* attribute), 42
- alarm\_77 (*AlarmsOverview* attribute), 46
- ALARM\_78 (*Alarms* attribute), 42
- alarm\_78 (*AlarmsOverview* attribute), 46
- ALARM\_79 (*Alarms* attribute), 42
- alarm\_79 (*AlarmsOverview* attribute), 46
- ALARM\_8 (*Alarms* attribute), 42
- alarm\_8 (*AlarmsOverview* attribute), 46
- ALARM\_80 (*Alarms* attribute), 42
- alarm\_80 (*AlarmsOverview* attribute), 46
- ALARM\_81 (*Alarms* attribute), 42
- alarm\_81 (*AlarmsOverview* attribute), 46
- ALARM\_82 (*Alarms* attribute), 42
- alarm\_82 (*AlarmsOverview* attribute), 46
- ALARM\_83 (*Alarms* attribute), 42
- alarm\_83 (*AlarmsOverview* attribute), 46
- ALARM\_84 (*Alarms* attribute), 42
- alarm\_84 (*AlarmsOverview* attribute), 46
- ALARM\_85 (*Alarms* attribute), 42
- alarm\_85 (*AlarmsOverview* attribute), 46
- ALARM\_86 (*Alarms* attribute), 42
- alarm\_86 (*AlarmsOverview* attribute), 46
- ALARM\_87 (*Alarms* attribute), 42
- alarm\_87 (*AlarmsOverview* attribute), 46
- ALARM\_88 (*Alarms* attribute), 42
- alarm\_88 (*AlarmsOverview* attribute), 46
- ALARM\_89 (*Alarms* attribute), 42
- alarm\_89 (*AlarmsOverview* attribute), 46
- ALARM\_9 (*Alarms* attribute), 42
- alarm\_9 (*AlarmsOverview* attribute), 46
- ALARM\_90 (*Alarms* attribute), 42
- alarm\_90 (*AlarmsOverview* attribute), 46
- ALARM\_91 (*Alarms* attribute), 42
- alarm\_91 (*AlarmsOverview* attribute), 46
- ALARM\_92 (*Alarms* attribute), 42
- alarm\_92 (*AlarmsOverview* attribute), 46
- ALARM\_93 (*Alarms* attribute), 42
- alarm\_93 (*AlarmsOverview* attribute), 46
- ALARM\_94 (*Alarms* attribute), 42
- alarm\_94 (*AlarmsOverview* attribute), 46
- ALARM\_95 (*Alarms* attribute), 42
- alarm\_95 (*AlarmsOverview* attribute), 46
- ALARM\_96 (*Alarms* attribute), 42
- alarm\_96 (*AlarmsOverview* attribute), 46
- ALARM\_97 (*Alarms* attribute), 42
- alarm\_97 (*AlarmsOverview* attribute), 46
- ALARM\_98 (*Alarms* attribute), 42
- alarm\_98 (*AlarmsOverview* attribute), 46
- ALARM\_99 (*Alarms* attribute), 42
- alarm\_99 (*AlarmsOverview* attribute), 46
- Alarms (*class* in *hvl\_ccb.dev.cube.constants*), 38
- AlarmsOverview (*class* in *hvl\_ccb.dev.cube.constants*), 42
- AlarmText (*class* in *hvl\_ccb.dev.supercube2015.constants*), 64
- ALL (*TiePieOscilloscopeAutoResolutionModes* attribute), 89
- all\_measurements (*LuminosMeasurementType* attribute), 177
- all\_measurements\_types() (*LuminosMeasurementType* class method), 177
- amplitude (*TiePieGeneratorConfig* property), 85
- ANALOG (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 135
- analog\_control (*FuGProbusVDIRegisters* property), 115
- ANY (*LabJack.DeviceType* attribute), 127
- ANY (*LJMCommunicationConfig.ConnectionType* attribute), 16
- ANY (*LJMCommunicationConfig.DeviceType* attribute), 16
- ANY (*TiePieOscilloscopeTriggerKind* attribute), 84
- ARBITRARY (*TiePieGeneratorSignalType* attribute), 86
- AsyncCommunicationProtocol (*class* in *hvl\_ccb.comm.base*), 11
- AsyncCommunicationProtocolConfig (*class* in *hvl\_ccb.comm.base*), 13
- ATM (*Pressure* attribute), 184
- ATMOSPHERE (*Pressure* attribute), 184
- attenuation (*CryLasAttenuator* property), 95
- AUTO (*EarthingStickOperatingStatus* attribute), 47
- AUTO (*LaudaProRp245eConfig.OperationModeEnum* attribute), 136
- AUTO (*RTO1024.TriggerModes* attribute), 162
- auto\_install\_mode (*T560* property), 56
- auto\_install\_mode (*T560Config* attribute), 57
- auto\_laser\_on (*CryLasLaserConfig* attribute), 100
- auto\_resolution\_mode (*TiePieOscilloscopeConfig* property), 89
- AutoInstallMode (*class* in *hvl\_ccb.dev.highland\_t560.base*), 53
- AutoNumberNameEnum (*class* in *hvl\_ccb.utils.enum*), 185

## B

B (*SupercubeOpcEndpoint* attribute), 71  
 BA (*NewportConfigCommands* attribute), 142  
 backlash\_compensation (*NewportSMC100PPConfig* attribute), 150  
 backup\_waveform() (*RTO1024* method), 162  
 bar (*PfeifferTPG.PressureUnits* attribute), 156  
 BAR (*Pressure* attribute), 184  
 barometric\_pressure (*LuminosMeasurementType* attribute), 177  
 base\_velocity (*NewportSMC100PPConfig* attribute), 150  
 BaseCube (class in *hvl\_ccb.dev.cube.base*), 33  
 BaseCubeConfiguration (class in *hvl\_ccb.dev.cube.base*), 36  
 BaseCubeOpcUaCommunication (class in *hvl\_ccb.dev.cube.base*), 37  
 BaseCubeOpcUaCommunicationConfig (class in *hvl\_ccb.dev.cube.base*), 37  
 BATH\_TEMP (*LaudaProRp245eCommand* attribute), 134  
 baudrate (*CryLasAttenuatorSerialCommunicationConfig* attribute), 96  
 baudrate (*CryLasLaserSerialCommunicationConfig* attribute), 102  
 baudrate (*FuGSerialCommunicationConfig* attribute), 118  
 baudrate (*HeinzingerSerialCommunicationConfig* attribute), 124  
 baudrate (*LuminosSerialCommunicationConfig* attribute), 178  
 baudrate (*MBW973SerialCommunicationConfig* attribute), 141  
 baudrate (*NewportSMC100PPSerialCommunicationConfig* attribute), 153  
 baudrate (*PfeifferTPGSerialCommunicationConfig* attribute), 160  
 baudrate (*SerialCommunicationConfig* attribute), 24  
 BH (*NewportConfigCommands* attribute), 142  
 board (*VisaCommunicationConfig* attribute), 31  
 breakdown\_detection\_active (*BaseCube* property), 33  
 breakdown\_detection\_reset() (*BaseCube* method), 34  
 breakdown\_detection\_triggered (*BaseCube* property), 34  
 BreakdownDetection (class in *hvl\_ccb.dev.supercube2015.constants*), 65  
 bufsize (*TcpCommunicationConfig* attribute), 27  
 build\_str() (*LaudaProRp245eCommand* method), 135  
 bytesize (*CryLasAttenuatorSerialCommunicationConfig* attribute), 96  
 bytesize (*CryLasLaserSerialCommunicationConfig* attribute), 103

bytesize (*FuGSerialCommunicationConfig* attribute), 118  
 bytesize (*HeinzingerSerialCommunicationConfig* attribute), 124  
 bytesize (*LuminosSerialCommunicationConfig* attribute), 178  
 bytesize (*MBW973SerialCommunicationConfig* attribute), 141  
 bytesize (*NewportSMC100PPSerialCommunicationConfig* attribute), 154  
 bytesize (*PfeifferTPGSerialCommunicationConfig* attribute), 160  
 Bytesize (*SerialCommunicationConfig* attribute), 24  
 bytesize (*SerialCommunicationConfig* attribute), 24

## C

C (*LabJack.TemperatureUnit* attribute), 127  
 C (*LabJack.ThermocoupleType* attribute), 128  
 C (*Temperature* attribute), 184  
 calibration\_factor (*CryLasLaserConfig* attribute), 100  
 calibration\_factor (*LEM4000S* attribute), 183  
 calibration\_mode (*FuGProbusVDIRegisters* property), 115  
 cc\_mode (*FuGProbusVDIRegisters* property), 115  
 CCException, 189  
 cee16 (*GeneralSockets* attribute), 67  
 cee16\_socket (*BaseCube* property), 34  
 CELSIUS (*Temperature* attribute), 184  
 ch\_a (*T560* property), 56  
 ch\_b (*T560* property), 56  
 ch\_c (*T560* property), 56  
 ch\_d (*T560* property), 56  
 channels\_enabled (*TiePieOscilloscope* property), 88  
 check\_for\_error() (*NewportSMC100PPSerialCommunication* method), 152  
 check\_generic\_type() (in module *hvl\_ccb.utils.typing*), 186  
 check\_master\_slave\_config() (*PSI9000* method), 104  
 chunk\_size (*VisaCommunicationConfig* attribute), 31  
 clean\_amplitude() (*TiePieGeneratorConfig* method), 85  
 clean\_auto\_resolution\_mode() (*TiePieOscilloscopeConfig* static method), 89  
 clean\_coupling() (*TiePieOscilloscopeChannelConfig* static method), 82  
 clean\_enabled() (*TiePieGeneratorConfig* static method), 85  
 clean\_enabled() (*TiePieOscilloscopeChannelConfig* static method), 82  
 clean\_frequency() (*TiePieGeneratorConfig* method), 85

- `clean_input_range()` (*TiePieOscilloscopeChannelConfig method*), 82
- `clean_offset()` (*TiePieGeneratorConfig method*), 85
- `clean_pre_sample_ratio()` (*TiePieOscilloscopeConfig method*), 90
- `clean_probe_offset()` (*TiePieOscilloscopeChannelConfig method*), 82
- `clean_record_length()` (*TiePieOscilloscopeConfig method*), 90
- `clean_resolution()` (*TiePieOscilloscopeConfig static method*), 90
- `clean_sample_frequency()` (*TiePieOscilloscopeConfig method*), 90
- `clean_signal_type()` (*TiePieGeneratorConfig static method*), 85
- `clean_trigger_enabled()` (*TiePieOscilloscopeChannelConfig static method*), 82
- `clean_trigger_hysteresis()` (*TiePieOscilloscopeChannelConfig method*), 82
- `clean_trigger_kind()` (*TiePieOscilloscopeChannelConfig static method*), 82
- `clean_trigger_level()` (*TiePieOscilloscopeChannelConfig method*), 82
- `clean_trigger_level_mode()` (*TiePieOscilloscopeChannelConfig static method*), 82
- `clean_trigger_timeout()` (*TiePieOscilloscopeConfig method*), 90
- `clean_values()` (*AsyncCommunicationProtocolConfig method*), 13
- `clean_values()` (*BaseCubeConfiguration method*), 36
- `clean_values()` (*CryLasAttenuatorConfig method*), 95
- `clean_values()` (*CryLasLaserConfig method*), 100
- `clean_values()` (*EmptyConfig method*), 93, 188
- `clean_values()` (*FuGConfig method*), 110
- `clean_values()` (*HeinzingerConfig method*), 120
- `clean_values()` (*ILS2TConfig method*), 172
- `clean_values()` (*LaudaProRp245eConfig method*), 136
- `clean_values()` (*LaudaProRp245eTcpCommunicationConfig method*), 137
- `clean_values()` (*LJMCommunicationConfig method*), 17
- `clean_values()` (*LuminoxConfig method*), 176
- `clean_values()` (*MBW973Config method*), 139
- `clean_values()` (*ModbusTcpCommunicationConfig method*), 19
- `clean_values()` (*NewportSMC100PPConfig method*), 150
- `clean_values()` (*OpcUaCommunicationConfig method*), 21
- `clean_values()` (*PfeifferTPGConfig method*), 158
- `clean_values()` (*PICubeConfiguration method*), 51
- `clean_values()` (*PSI9000Config method*), 106
- `clean_values()` (*SerialCommunicationConfig method*), 24
- `clean_values()` (*SupercubeConfiguration method*), 62
- `clean_values()` (*T560Config method*), 57
- `clean_values()` (*TcpCommunicationConfig method*), 27
- `clean_values()` (*TechnixConfig method*), 78
- `clean_values()` (*TelnetCommunicationConfig method*), 29
- `clean_values()` (*TiePieDeviceConfig method*), 80
- `clean_values()` (*VisaCommunicationConfig method*), 31
- `clean_waveform()` (*TiePieGeneratorConfig method*), 85
- `Client` (*class in hvl\_ccb.comm.opc*), 20
- `CLOSE` (*EarthingStickOperation attribute*), 47
- `close()` (*CommunicationProtocol method*), 14
- `close()` (*LaudaProRp245eTcpCommunication method*), 137
- `close()` (*LJMCommunication method*), 15
- `close()` (*ModbusTcpCommunication method*), 18
- `close()` (*NullCommunicationProtocol method*), 14
- `close()` (*OpcUaCommunication method*), 20
- `close()` (*SerialCommunication method*), 23
- `close()` (*Tcp method*), 26
- `close()` (*TelnetCommunication method*), 28
- `close()` (*VisaCommunication method*), 30
- `close_shutter()` (*CryLasLaser method*), 98
- `CLOSED` (*CryLasLaser.AnswersShutter attribute*), 97
- `CLOSED` (*CryLasLaserShutterStatus attribute*), 103
- `CLOSED` (*DoorStatus attribute*), 47
- `closed` (*DoorStatus attribute*), 65
- `CLOSED` (*EarthingStickStatus attribute*), 48
- `closed` (*EarthingStickStatus attribute*), 66
- `CMD_EXEC_ERROR` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- `CMD_NOT_ALLOWED` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- `CMD_NOT_ALLOWED_CC` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- `CMD_NOT_ALLOWED_CONFIGURATION` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- `CMD_NOT_ALLOWED_DISABLE` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- `CMD_NOT_ALLOWED_HOMING` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- `CMD_NOT_ALLOWED_MOVING` (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151

attribute), 151

CMD\_NOT\_ALLOWED\_NOT\_REFERENCED (New-  
portSMC100PPSerialCommunication.ControllerErrors  
attribute), 151

CMD\_NOT\_ALLOWED\_PP (New-  
portSMC100PPSerialCommunication.ControllerErrors  
attribute), 151

CMD\_NOT\_ALLOWED\_READY (New-  
portSMC100PPSerialCommunication.ControllerErrors  
attribute), 151

CMR (PfeifferTPG.SensorTypes attribute), 156

CODE\_OR\_ADDR\_INVALID (New-  
portSMC100PPSerialCommunication.ControllerErrors  
attribute), 151

collect\_measurement\_data() (TiePieOscilloscope  
method), 88

com (SingleCommDevice property), 94

COM\_TIME\_OUT (LaudaProRp245eCommand attribute),  
134

com\_time\_out (LaudaProRp245eConfig attribute), 136

COM\_TIMEOUT (NewportSMC100PPSerialCommunication.ControllerErrors  
attribute), 151

command (LuminoxMeasurementType property), 177

COMMAND (TriggerMode attribute), 55

command() (FuGProbusIV method), 113

communication\_channel (TechnixConfig attribute), 78

CommunicationException, 14

CommunicationProtocol (class in  
hvl\_ccb.comm.base), 14

config (ConfigurationMixin property), 187

CONFIG (FuGProbusVRegisterGroups attribute), 116

CONFIG (NewportSMC100PP.StateMessages attribute),  
143

CONFIG (NewportStates attribute), 154

config\_cls() (AsyncCommunicationProtocol static  
method), 11

config\_cls() (BaseCube static method), 34

config\_cls() (BaseCubeOpcUaCommunication static  
method), 37

config\_cls() (ConfigurationMixin static method), 187

config\_cls() (CryLasAttenuator static method), 95

config\_cls() (CryLasAttenuatorSerialCommunication  
static method), 96

config\_cls() (CryLasLaser static method), 98

config\_cls() (CryLasLaserSerialCommunication  
static method), 102

config\_cls() (Device static method), 92

config\_cls() (FuGProbusIV static method), 113

config\_cls() (FuGSerialCommunication static  
method), 118

config\_cls() (HeinzingerDI static method), 121

config\_cls() (HeinzingerSerialCommunication static  
method), 124

config\_cls() (ILS2T static method), 170

config\_cls() (ILS2TModbusTcpCommunication static  
method), 173

config\_cls() (LaudaProRp245e static method), 132

config\_cls() (LaudaProRp245eTcpCommunication  
static method), 137

config\_cls() (LJMCommunication static method), 15

config\_cls() (Luminox static method), 175

config\_cls() (LuminoxSerialCommunication static  
method), 178

config\_cls() (MBW973 static method), 138

config\_cls() (MBW973SerialCommunication static  
method), 140

config\_cls() (ModbusTcpCommunication static  
method), 18

config\_cls() (NewportSMC100PP static method), 144

config\_cls() (NewportSMC100PPSerialCommunication  
static method), 152

config\_cls() (NullCommunicationProtocol static  
method), 14

config\_cls() (OpcUaCommunication static method),  
159

config\_cls() (PfeifferTPG static method), 156

config\_cls() (PfeifferTPGSerialCommunication static  
method), 159

config\_cls() (PICube static method), 50

config\_cls() (PICubeOpcUaCommunication static  
method), 51

config\_cls() (PSI9000 static method), 104

config\_cls() (PSI9000VisaCommunication static  
method), 107

config\_cls() (RTO1024 static method), 162

config\_cls() (RTO1024VisaCommunication static  
method), 168

config\_cls() (SerialCommunication static method), 23

config\_cls() (Supercube2015Base static method), 58

config\_cls() (SupercubeAOPcUaCommunication  
static method), 72

config\_cls() (SupercubeOpcUaCommunication static  
method), 62

config\_cls() (SyncCommunicationProtocol static  
method), 14

config\_cls() (T560 static method), 56

config\_cls() (T560Communication static method), 54

config\_cls() (Tcp static method), 26

config\_cls() (Technix static method), 77

config\_cls() (TechnixSerialCommunication static  
method), 74

config\_cls() (TechnixTelnetCommunication static  
method), 75

config\_cls() (TelnetCommunication static method), 28

config\_cls() (TiePieOscilloscope static method), 88

config\_cls() (VisaCommunication static method), 30

config\_cls() (VisaDevice static method), 181

config\_gen (TiePieGeneratorMixin attribute), 86

- config\_high\_pulse() (*LabJack method*), 128  
 config\_i2c (*TiePieI2CHostMixin attribute*), 87  
 config\_osc (*TiePieOscilloscope attribute*), 88  
 config\_osc\_channel\_dict (*TiePieOscilloscope attribute*), 88  
 config\_status (*FuG property*), 108  
 configdataclass() (*in module hvl\_ccb.configuration*), 188  
 configuration\_save\_json() (*ConfigurationMixin method*), 187  
 ConfigurationMixin (*class in hvl\_ccb.configuration*), 187  
 ConfigurationValueWarning, 187  
 connection\_type (*LJMCommunicationConfig attribute*), 17  
 CONT\_MODE (*LaudaProRp245eCommand attribute*), 134  
 continue\_ramp() (*LaudaProRp245e method*), 132  
 control\_mode (*LaudaProRp245eConfig attribute*), 136  
 CONVERSION (*LEM4000S attribute*), 182  
 convert() (*LEM4000S method*), 183  
 convert() (*LMT70A method*), 183  
 convert() (*Sensor method*), 183  
 convert() (*Unit class method*), 184  
 ConvertableTypes (*in module hvl\_ccb.utils.typing*), 186  
 COOLOFF (*LaudaProRp245eConfig.OperationModeEnum attribute*), 136  
 COOLON (*LaudaProRp245eConfig.OperationModeEnum attribute*), 136  
 coupling (*TiePieOscilloscopeChannelConfig property*), 82  
 CR (*FuGTerminators attribute*), 119  
 create\_serial\_port() (*SerialCommunicationConfig method*), 24  
 create\_telnet() (*TelnetCommunicationConfig method*), 29  
 CRLF (*FuGTerminators attribute*), 119  
 CryLasAttenuator (*class in hvl\_ccb.dev.crylas*), 95  
 CryLasAttenuatorConfig (*class in hvl\_ccb.dev.crylas*), 95  
 CryLasAttenuatorError, 96  
 CryLasAttenuatorSerialCommunication (*class in hvl\_ccb.dev.crylas*), 96  
 CryLasAttenuatorSerialCommunicationConfig (*class in hvl\_ccb.dev.crylas*), 96  
 CryLasLaser (*class in hvl\_ccb.dev.crylas*), 97  
 CryLasLaser.AnswersShutter (*class in hvl\_ccb.dev.crylas*), 97  
 CryLasLaser.AnswersStatus (*class in hvl\_ccb.dev.crylas*), 97  
 CryLasLaser.LaserStatus (*class in hvl\_ccb.dev.crylas*), 98  
 CryLasLaser.RepetitionRates (*class in hvl\_ccb.dev.crylas*), 98  
 CryLasLaserConfig (*class in hvl\_ccb.dev.crylas*), 100  
 CryLasLaserError, 101  
 CryLasLaserNotReadyError, 101  
 CryLasLaserPoller (*class in hvl\_ccb.dev.crylas*), 101  
 CryLasLaserSerialCommunication (*class in hvl\_ccb.dev.crylas*), 101  
 CryLasLaserSerialCommunicationConfig (*class in hvl\_ccb.dev.crylas*), 102  
 CryLasLaserShutterStatus (*class in hvl\_ccb.dev.crylas*), 103  
 CubeEarthingStickOperationError, 46  
 CubeError, 47  
 CubeRemoteControlError, 47  
 CubeStatusChangeError, 47  
 CubeStopError, 47  
 current (*FuG property*), 108  
 CURRENT (*FuGProbusIVCommands attribute*), 113  
 CURRENT (*FuGReadbackChannels attribute*), 117  
 current (*Technix property*), 77  
 current\_lower\_limit (*PSI9000Config attribute*), 106  
 current\_monitor (*FuG property*), 109  
 current\_primary (*PICube property*), 50  
 current\_primary (*Power attribute*), 69  
 current\_upper\_limit (*PSI9000Config attribute*), 106  
 CurrentPosition (*New-portSMC100PPConfig.HomeSearch attribute*), 149  
 cv\_mode (*FuGProbusVDIRegisters property*), 115
- ## D
- datachange\_notification() (*OpcUaSubHandler method*), 22  
 datachange\_notification() (*SupercubeSubscriptionHandler method*), 63  
 date\_of\_manufacture (*LuminosMeasurementType attribute*), 177  
 datetime\_to\_opc() (*BaseCube class method*), 34  
 DC (*TiePieGeneratorSignalType attribute*), 86  
 DC\_140KV (*PowerSetup attribute*), 49  
 DC\_280KV (*PowerSetup attribute*), 49  
 DC\_DoubleStage\_280kV (*PowerSetup attribute*), 70  
 DC\_SingleStage\_140kV (*PowerSetup attribute*), 70  
 DC\_VOLTAGE\_TOO\_LOW (*New-portSMC100PP.MotorErrors attribute*), 143  
 DCA (*TiePieOscilloscopeChannelCoupling attribute*), 83  
 DCV (*TiePieOscilloscopeChannelCoupling attribute*), 83  
 default\_com\_cls() (*BaseCube static method*), 34  
 default\_com\_cls() (*CryLasAttenuator static method*), 95  
 default\_com\_cls() (*CryLasLaser static method*), 98  
 default\_com\_cls() (*FuGProbusIV static method*), 113  
 default\_com\_cls() (*HeinzingerDI static method*), 121  
 default\_com\_cls() (*ILS2T static method*), 170

- default\_com\_cls() (*LabJack static method*), 128
- default\_com\_cls() (*LaudaProRp245e static method*), 132
- default\_com\_cls() (*Luminox static method*), 176
- default\_com\_cls() (*MBW973 static method*), 138
- default\_com\_cls() (*NewportSMC100PP static method*), 144
- default\_com\_cls() (*PfeifferTPG static method*), 156
- default\_com\_cls() (*PICube static method*), 50
- default\_com\_cls() (*PSI9000 static method*), 104
- default\_com\_cls() (*RTO1024 static method*), 162
- default\_com\_cls() (*SingleCommDevice static method*), 94
- default\_com\_cls() (*Supercube2015Base static method*), 59
- default\_com\_cls() (*Supercube2015WithFU static method*), 71
- default\_com\_cls() (*T560 static method*), 56
- default\_com\_cls() (*Technix method*), 77
- default\_com\_cls() (*TiePieOscilloscope static method*), 88
- default\_com\_cls() (*VisaDevice static method*), 181
- DEFAULT\_IO\_SCANNING\_CONTROL\_VALUES (*ILS2T attribute*), 169
- default\_n\_attempts\_read\_text\_nonempty (*AsyncCommunicationProtocolConfig attribute*), 13
- default\_n\_attempts\_read\_text\_nonempty (*FuGSerialCommunicationConfig attribute*), 118
- default\_n\_attempts\_read\_text\_nonempty (*HeinzingerSerialCommunicationConfig attribute*), 124
- default\_number\_of\_recordings (*HeinzingerConfig attribute*), 120
- Device (*class in hvl\_ccb.dev.base*), 92
- DEVICE\_TYPE (*LaudaProRp245eCommand attribute*), 134
- device\_type (*LJMCommunicationConfig attribute*), 17
- DeviceException, 92
- DeviceExistingException, 92
- DeviceFailuresException, 92
- devices\_failed\_start (*DeviceSequenceMixin attribute*), 93
- devices\_failed\_start (*ExperimentManager attribute*), 190
- devices\_failed\_stop (*DeviceSequenceMixin attribute*), 93
- devices\_failed\_stop (*ExperimentManager attribute*), 190
- DeviceSequenceMixin (*class in hvl\_ccb.dev.base*), 92
- di (*FuG property*), 109
- digital\_control (*FuGProbusVDIRegisters property*), 115
- DIOChannel (*LabJack attribute*), 127
- DISABLE (*NewportStates attribute*), 154
- disable() (*ILS2T method*), 170
- DISABLE\_FROM\_JOGGING (*NewportSMC100PP.StateMessages attribute*), 143
- DISABLE\_FROM\_MOVING (*NewportSMC100PP.StateMessages attribute*), 143
- DISABLE\_FROM\_READY (*NewportSMC100PP.StateMessages attribute*), 143
- disable\_pulses() (*LabJack method*), 128
- DISABLED (*TiePieOscilloscopeAutoResolutionModes attribute*), 89
- DisableEspStageCheck (*NewportSMC100PPConfig.EspStageConfig attribute*), 149
- disarm\_trigger() (*T560 method*), 57
- disconnect() (*Client method*), 20
- DISPLACEMENT\_OUT\_OF\_LIMIT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 151
- display\_message\_board() (*BaseCube method*), 34
- display\_status\_board() (*BaseCube method*), 34
- do\_ioscanning\_write() (*ILS2T method*), 170
- door\_1\_status (*BaseCube attribute*), 34
- door\_2\_status (*BaseCube attribute*), 34
- door\_3\_status (*BaseCube attribute*), 34
- DoorStatus (*class in hvl\_ccb.dev.cube.constants*), 47
- DoorStatus (*class in hvl\_ccb.dev.supercube2015.constants*), 65

## E

- E (*LabJack.ThermocoupleType attribute*), 128
- E0 (*FuGErrorcodes attribute*), 111
- E1 (*FuGErrorcodes attribute*), 111
- E10 (*FuGErrorcodes attribute*), 111
- E100 (*FuGErrorcodes attribute*), 111
- E106 (*FuGErrorcodes attribute*), 111
- E11 (*FuGErrorcodes attribute*), 111
- E115 (*FuGErrorcodes attribute*), 111
- E12 (*FuGErrorcodes attribute*), 111
- E125 (*FuGErrorcodes attribute*), 111
- E13 (*FuGErrorcodes attribute*), 111
- E135 (*FuGErrorcodes attribute*), 111
- E14 (*FuGErrorcodes attribute*), 111
- E145 (*FuGErrorcodes attribute*), 111
- E15 (*FuGErrorcodes attribute*), 111
- E155 (*FuGErrorcodes attribute*), 111
- E16 (*FuGErrorcodes attribute*), 112
- E165 (*FuGErrorcodes attribute*), 112
- E2 (*FuGErrorcodes attribute*), 112
- E206 (*FuGErrorcodes attribute*), 112
- E306 (*FuGErrorcodes attribute*), 112
- E4 (*FuGErrorcodes attribute*), 112

E5 (*FuErrorcodes* attribute), 112  
E504 (*FuErrorcodes* attribute), 112  
E505 (*FuErrorcodes* attribute), 112  
E6 (*FuErrorcodes* attribute), 112  
E666 (*FuErrorcodes* attribute), 112  
E7 (*FuErrorcodes* attribute), 112  
E8 (*FuErrorcodes* attribute), 112  
E9 (*FuErrorcodes* attribute), 112  
earthing\_rod\_1\_status (*BaseCube* attribute), 34  
earthing\_rod\_2\_status (*BaseCube* attribute), 34  
earthing\_rod\_3\_status (*BaseCube* attribute), 34  
EarthingRodStatus (class *hvl\_ccb.dev.cube.constants*), 47  
EarthingStick (class *hvl\_ccb.dev.supercube2015.constants*), 65  
EarthingStickOperatingStatus (class *hvl\_ccb.dev.cube.constants*), 47  
EarthingStickOperation (class *hvl\_ccb.dev.cube.constants*), 47  
EarthingStickStatus (class *hvl\_ccb.dev.cube.constants*), 48  
EarthingStickStatus (class *hvl\_ccb.dev.supercube2015.constants*), 66  
EEPROM\_ACCESS\_ERROR (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 151  
EIGHT (*HeinzingerConfig.RecordingsEnum* attribute), 120  
EIGHT\_BIT (*TiePieOscilloscopeResolution* attribute), 90  
EIGHT\_HUNDRED\_MILLI\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
EIGHT\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
EIGHTBITS (*SerialCommunicationBytesize* attribute), 24  
EIGHTY\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
EmptyConfig (class in *hvl\_ccb.configuration*), 188  
EmptyConfig (class in *hvl\_ccb.dev.base*), 93  
enable() (*ILS2T* method), 170  
enable\_clock() (*LabJack* method), 128  
enabled (*TiePieGeneratorConfig* property), 85  
enabled (*TiePieOscilloscopeChannelConfig* property), 82  
EnableEspStageCheck (*NewportSMC100PPConfig.EspStageConfig* attribute), 149  
encoding (*AsyncCommunicationProtocolConfig* attribute), 13  
encoding (*NewportSMC100PPSerialCommunicationConfig* attribute), 154  
encoding\_error\_handling (*AsyncCommunicationProtocolConfig* attribute), 13  
encoding\_error\_handling (*NewportSMC100PPSerialCommunicationConfig* attribute), 154  
EndOfRunSwitch (*NewportSMC100PPConfig.HomeSearch* attribute), 149  
EndOfRunSwitch\_and\_Index (*NewportSMC100PPConfig.HomeSearch* attribute), 149  
endpoint\_name (*BaseCubeOpcUaCommunicationConfig* attribute), 37  
endpoint\_name (*OpcUaCommunicationConfig* attribute), 21  
endpoint\_name (*PICubeOpcUaCommunicationConfig* attribute), 52  
in endpoint\_name (*SupercubeAOpcUaConfiguration* attribute), 73  
in ERROR (*DoorStatus* attribute), 47  
error (*DoorStatus* attribute), 65  
in ERROR (*EarthingStickStatus* attribute), 48  
error (*EarthingStickStatus* attribute), 67  
in ERROR (*ExperimentStatus* attribute), 190  
ERROR (*SafetyStatus* attribute), 49  
in Error (*SafetyStatus* attribute), 70  
errorcode (*FuGError* attribute), 111  
in Errors (class in *hvl\_ccb.dev.supercube2015.constants*), 67  
ESP\_STAGE\_NAME\_INVALID (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 151  
ETH (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 135  
ETHERNET (*LJMCommunicationConfig.ConnectionType* attribute), 16  
EVEN (*SerialCommunicationParity* attribute), 25  
event\_notification() (*OpcUaSubHandler* method), 22  
EXECUTE (*FuGProbusIVCommands* attribute), 113  
execute\_absolute\_position() (*ILS2T* method), 170  
execute\_on\_x (*FuGProbusVConfigRegisters* property), 114  
execute\_relative\_step() (*ILS2T* method), 170  
EXECUTEONX (*FuGProbusIVCommands* attribute), 113  
exit\_configuration() (*NewportSMC100PP* method), 144  
exit\_configuration\_wait\_sec (*NewportSMC100PPConfig* attribute), 150  
EXPERIMENT\_BLOCKED (*EarthingRodStatus* attribute), 47  
EXPERIMENT\_READY (*EarthingRodStatus* attribute), 47  
ExperimentError, 189  
ExperimentManager (class in *hvl\_ccb.experiment\_manager*), 189  
ExperimentStatus (class in *hvl\_ccb.experiment\_manager*), 190  
EXPT100 (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 135  
EXT\_FALLING\_EDGE (*TriggerMode* attribute), 55  
EXT\_RISING\_EDGE (*TriggerMode* attribute), 55

- External (*PowerSetup* attribute), 70  
 EXTERNAL\_SOURCE (*PowerSetup* attribute), 49  
 EXTERNAL\_TEMP (*LaudaProRp245eCommand* attribute), 134
- ## F
- F (*LabJack.TemperatureUnit* attribute), 127  
 F (*Temperature* attribute), 184  
 FAHRENHEIT (*Temperature* attribute), 184  
 failures (*DeviceFailuresException* attribute), 92  
 FALLING (*TiePieOscilloscopeTriggerKind* attribute), 84  
 FAST (*ILS2T.Ref16Jog* attribute), 169  
 file\_copy() (*RTO1024* method), 162  
 finish() (*ExperimentManager* method), 190  
 FINISHED (*ExperimentStatus* attribute), 190  
 FINISHING (*ExperimentStatus* attribute), 190  
 fire\_trigger() (*T560* method), 57  
 FIRMWARE (*FuGReadbackChannels* attribute), 117  
 FIVE\_MHZ (*LabJack.ClockFrequency* attribute), 126  
 FIVEBITS (*SerialCommunicationBytesize* attribute), 24  
 FLT\_INFO (*ILS2TRegAddr* attribute), 174  
 FLT\_MEM\_DEL (*ILS2TRegAddr* attribute), 174  
 FLT\_MEM\_RESET (*ILS2TRegAddr* attribute), 174  
 FOLLOWING\_ERROR (*NewportSMC100PP.MotorErrors* attribute), 143  
 FOLLOWRAMP (*FuGRampModes* attribute), 117  
 force\_trigger() (*TiePieOscilloscope* method), 89  
 force\_value() (*AsyncCommunicationProtocolConfig* method), 13  
 force\_value() (*BaseCubeConfiguration* method), 36  
 force\_value() (*BaseCubeOpcUaCommunicationConfig* method), 37  
 force\_value() (*CryLasAttenuatorConfig* method), 95  
 force\_value() (*CryLasAttenuatorSerialCommunicationConfig* method), 96  
 force\_value() (*CryLasLaserConfig* method), 100  
 force\_value() (*CryLasLaserSerialCommunicationConfig* method), 103  
 force\_value() (*EmptyConfig* method), 93, 188  
 force\_value() (*FuGConfig* method), 110  
 force\_value() (*FuGSerialCommunicationConfig* method), 118  
 force\_value() (*HeinzingerConfig* method), 120  
 force\_value() (*HeinzingerSerialCommunicationConfig* method), 124  
 force\_value() (*ILS2TConfig* method), 172  
 force\_value() (*ILS2TModbusTcpCommunicationConfig* method), 173  
 force\_value() (*LaudaProRp245eConfig* method), 136  
 force\_value() (*LaudaProRp245eTcpCommunicationConfig* method), 137  
 force\_value() (*LJMCommunicationConfig* method), 17  
 force\_value() (*LuminoxConfig* method), 176  
 force\_value() (*LuminoxSerialCommunicationConfig* method), 179  
 force\_value() (*MBW973Config* method), 140  
 force\_value() (*MBW973SerialCommunicationConfig* method), 141  
 force\_value() (*ModbusTcpCommunicationConfig* method), 19  
 force\_value() (*NewportSMC100PPConfig* method), 150  
 force\_value() (*NewportSMC100PPSerialCommunicationConfig* method), 154  
 force\_value() (*OpcUaCommunicationConfig* method), 21  
 force\_value() (*PfeifferTPGConfig* method), 158  
 force\_value() (*PfeifferTPGSerialCommunicationConfig* method), 160  
 force\_value() (*PICubeConfiguration* method), 51  
 force\_value() (*PICubeOpcUaCommunicationConfig* method), 52  
 force\_value() (*PSI9000Config* method), 106  
 force\_value() (*PSI9000VisaCommunicationConfig* method), 107  
 force\_value() (*RTO1024Config* method), 167  
 force\_value() (*RTO1024VisaCommunicationConfig* method), 168  
 force\_value() (*SerialCommunicationConfig* method), 24  
 force\_value() (*SupercubeAOpcUaConfiguration* method), 73  
 force\_value() (*SupercubeConfiguration* method), 62  
 force\_value() (*SupercubeOpcUaCommunicationConfig* method), 63  
 force\_value() (*T560CommunicationConfig* method), 54  
 force\_value() (*T560Config* method), 57  
 force\_value() (*TcpCommunicationConfig* method), 27  
 force\_value() (*TechnixConfig* method), 78  
 force\_value() (*TechnixSerialCommunicationConfig* method), 75  
 force\_value() (*TechnixTelnetCommunicationConfig* method), 76  
 force\_value() (*TelnetCommunicationConfig* method), 29  
 force\_value() (*TiePieDeviceConfig* method), 80  
 force\_value() (*VisaCommunicationConfig* method), 31  
 force\_value() (*VisaDeviceConfig* method), 181  
 FORTY\_MHZ (*LabJack.ClockFrequency* attribute), 127  
 FORTY\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
 FOUR (*HeinzingerConfig.RecordingsEnum* attribute), 120  
 FOUR\_HUNDRED\_MILLI\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
 FOUR\_VOLT (*TiePieOscilloscopeRange* attribute), 83

FOURTEEN\_BIT (*TiePieOscilloscopeResolution* attribute), 90

FREERUN (*RTO1024.TriggerModes* attribute), 162

frequency (*PICube* property), 50

frequency (*Power* attribute), 69

frequency (*T560* property), 57

frequency (*TiePieGeneratorConfig* property), 85

FRM (*NewportConfigCommands* attribute), 142

from\_json() (*ConfigurationMixin* class method), 187

FRS (*NewportConfigCommands* attribute), 142

fso\_reset() (*Supercube2015WithFU* method), 72

FuG (class in *hvl\_ccb.dev.fug*), 108

FuGConfig (class in *hvl\_ccb.dev.fug*), 110

FuGDigitalVal (class in *hvl\_ccb.dev.fug*), 110

FuGError, 111

FuGErrorcodes (class in *hvl\_ccb.dev.fug*), 111

FuGMonitorModes (class in *hvl\_ccb.dev.fug*), 112

FuGPolarities (class in *hvl\_ccb.dev.fug*), 113

FuGProbusIV (class in *hvl\_ccb.dev.fug*), 113

FuGProbusIVCommands (class in *hvl\_ccb.dev.fug*), 113

FuGProbusV (class in *hvl\_ccb.dev.fug*), 114

FuGProbusVConfigRegisters (class in *hvl\_ccb.dev.fug*), 114

FuGProbusVDIRegisters (class in *hvl\_ccb.dev.fug*), 115

FuGProbusVDORegisters (class in *hvl\_ccb.dev.fug*), 115

FuGProbusVMonitorRegisters (class in *hvl\_ccb.dev.fug*), 116

FuGProbusVRegisterGroups (class in *hvl\_ccb.dev.fug*), 116

FuGProbusVSetRegisters (class in *hvl\_ccb.dev.fug*), 116

FuGRampModes (class in *hvl\_ccb.dev.fug*), 117

FuGReadbackChannels (class in *hvl\_ccb.dev.fug*), 117

FuGSerialCommunication (class in *hvl\_ccb.dev.fug*), 117

FuGSerialCommunicationConfig (class in *hvl\_ccb.dev.fug*), 118

FuGTerminators (class in *hvl\_ccb.dev.fug*), 119

**G**

gate\_mode (*T560* property), 57

gate\_polarity (*T560* property), 57

GateMode (class in *hvl\_ccb.dev.highland\_t560.base*), 53

GeneralSockets (class in *hvl\_ccb.dev.supercube2015.constants*), 67

GeneralSupport (class in *hvl\_ccb.dev.supercube2015.constants*), 67

GENERATOR (*TiePieDeviceType* attribute), 81

generator\_start() (*TiePieGeneratorMixin* method), 86

generator\_stop() (*TiePieGeneratorMixin* method), 86

get() (*AlarmText* class method), 64

get() (*MeasurementsDividerRatio* class method), 68

get() (*MeasurementsScaledInput* class method), 69

get\_acceleration() (*NewportSMC100PP* method), 144

get\_acquire\_length() (*RTO1024* method), 162

get\_ain() (*LabJack* method), 128

get\_bath\_temp() (*LaudaProRp245e* method), 132

get\_by\_p\_id() (*LabJack.DeviceType* class method), 127

get\_by\_p\_id() (*LJMCommunicationConfig.DeviceType* class method), 16

get\_cal\_current\_source() (*LabJack* method), 128

get\_cee16\_socket() (*Supercube2015Base* method), 59

get\_channel\_offset() (*RTO1024* method), 162

get\_channel\_position() (*RTO1024* method), 163

get\_channel\_range() (*RTO1024* method), 163

get\_channel\_scale() (*RTO1024* method), 163

get\_channel\_state() (*RTO1024* method), 163

get\_clock() (*LabJack* method), 129

get\_controller\_information() (*NewportSMC100PP* method), 144

get\_current() (*HeinzingerDI* method), 121

get\_dc\_volt() (*ILS2T* method), 171

get\_device() (*DeviceSequenceMixin* method), 93

get\_device\_by\_serial\_number() (in module *hvl\_ccb.dev.tiepie.base*), 81

get\_device\_type() (*LaudaProRp245e* method), 132

get\_devices() (*DeviceSequenceMixin* method), 93

get\_digital\_input() (*LabJack* method), 129

get\_door\_status() (*Supercube2015Base* method), 59

get\_earthing\_manual() (*Supercube2015Base* method), 59

get\_earthing\_status() (*Supercube2015Base* method), 59

get\_error\_code() (*ILS2T* method), 171

get\_error\_queue() (*VisaDevice* method), 181

get\_frequency() (*Supercube2015WithFU* method), 72

get\_fso\_active() (*Supercube2015WithFU* method), 72

get\_full\_scale\_mbar() (*PfeifferTPG* method), 156

get\_full\_scale\_unitless() (*PfeifferTPG* method), 157

get\_identification() (*VisaDevice* method), 181

get\_interface\_version() (*HeinzingerDI* method), 121

get\_max\_voltage() (*Supercube2015WithFU* method), 72

get\_measurement\_ratio() (*Supercube2015Base* method), 59

get\_measurement\_voltage() (*Supercube2015Base* method), 59

get\_motor\_configuration() (*NewportSMC100PP* method), 145

- get\_move\_duration() (*NewportSMC100PP method*), 145
- get\_negative\_software\_limit() (*NewportSMC100PP method*), 145
- get\_number\_of\_recordings() (*HeinzingerDI method*), 121
- get\_objects\_node() (*Client method*), 20
- get\_objects\_node() (*Server method*), 22
- get\_output() (*PSI9000 method*), 104
- get\_position() (*ILS2T method*), 171
- get\_position() (*NewportSMC100PP method*), 145
- get\_positive\_software\_limit() (*NewportSMC100PP method*), 145
- get\_power\_setup() (*Supercube2015WithFU method*), 72
- get\_primary\_current() (*Supercube2015WithFU method*), 72
- get\_primary\_voltage() (*Supercube2015WithFU method*), 72
- get\_product\_id() (*LabJack method*), 129
- get\_product\_name() (*LabJack method*), 129
- get\_product\_type() (*LabJack method*), 129
- get\_pulse\_energy\_and\_rate() (*CryLasLaser method*), 98
- get\_reference\_point() (*RTO1024 method*), 163
- get\_register() (*FuGProbusV method*), 114
- get\_repetitions() (*RTO1024 method*), 163
- get\_sbus\_rh() (*LabJack method*), 129
- get\_sbus\_temp() (*LabJack method*), 129
- get\_serial\_number() (*HeinzingerDI method*), 121
- get\_serial\_number() (*LabJack method*), 129
- get\_state() (*NewportSMC100PP method*), 146
- get\_status() (*ILS2T method*), 171
- get\_status() (*Supercube2015Base method*), 59
- get\_support\_input() (*Supercube2015Base method*), 59
- get\_support\_output() (*Supercube2015Base method*), 60
- get\_system\_lock() (*PSI9000 method*), 104
- get\_t13\_socket() (*Supercube2015Base method*), 60
- get\_target\_voltage() (*Supercube2015WithFU method*), 72
- get\_temperature() (*ILS2T method*), 171
- get\_timestamps() (*RTO1024 method*), 163
- get\_ui\_lower\_limits() (*PSI9000 method*), 105
- get\_uip\_upper\_limits() (*PSI9000 method*), 105
- get\_voltage() (*HeinzingerDI method*), 121
- get\_voltage\_current\_setpoint() (*PSI9000 method*), 105
- go\_home() (*NewportSMC100PP method*), 146
- go\_to\_configuration() (*NewportSMC100PP method*), 146
- GREEN\_NOT\_READY (*SafetyStatus attribute*), 49
- GREEN\_READY (*SafetyStatus attribute*), 49
- GreenNotReady (*SafetyStatus attribute*), 70
- GreenReady (*SafetyStatus attribute*), 70
- ## H
- HARDWARE (*CryLasLaser.RepetitionRates attribute*), 98
- has\_safe\_ground (*TiePieOscilloscopeChannelConfig property*), 82
- HEAD (*CryLasLaser.AnswersStatus attribute*), 97
- HeinzingerConfig (*class in hvl\_ccb.dev.heinzinger*), 119
- HeinzingerConfig.RecordingsEnum (*class in hvl\_ccb.dev.heinzinger*), 120
- HeinzingerDI (*class in hvl\_ccb.dev.heinzinger*), 120
- HeinzingerDI.OutputStatus (*class in hvl\_ccb.dev.heinzinger*), 121
- HeinzingerPNC (*class in hvl\_ccb.dev.heinzinger*), 122
- HeinzingerPNC.UnitCurrent (*class in hvl\_ccb.dev.heinzinger*), 122
- HeinzingerPNC.UnitVoltage (*class in hvl\_ccb.dev.heinzinger*), 122
- HeinzingerPNCDeviceNotRecognizedException, 123
- HeinzingerPNCError, 123
- HeinzingerPNCMaxCurrentExceededException, 123
- HeinzingerPNCMaxVoltageExceededException, 123
- HeinzingerSerialCommunication (*class in hvl\_ccb.dev.heinzinger*), 123
- HeinzingerSerialCommunicationConfig (*class in hvl\_ccb.dev.heinzinger*), 124
- HIGH (*LabJack.DIOStatus attribute*), 127
- high\_resolution (*FuGProbusVSetRegisters property*), 116
- home\_search\_polling\_interval (*NewportSMC100PPConfig attribute*), 150
- home\_search\_timeout (*NewportSMC100PPConfig attribute*), 150
- home\_search\_type (*NewportSMC100PPConfig attribute*), 150
- home\_search\_velocity (*NewportSMC100PPConfig attribute*), 150
- HOME\_STARTED (*NewportSMC100PPSerialCommunication.ControllerError attribute*), 151
- HomeSwitch (*NewportSMC100PPConfig.HomeSearch attribute*), 149
- HomeSwitch\_and\_Index (*NewportSMC100PPConfig.HomeSearch attribute*), 149
- HOMING (*NewportStates attribute*), 154
- HOMING\_FROM\_RS232 (*NewportSMC100PP.StateMessages attribute*), 143
- HOMING\_FROM\_SMC (*NewportSMC100PP.StateMessages attribute*), 143

HOMING\_TIMEOUT (*NewportSMC100PP.MotorErrors attribute*), 143  
 horn (*Safety attribute*), 70  
 horn() (*Supercube2015Base method*), 60  
 host (*ModbusTcpCommunicationConfig attribute*), 19  
 host (*OpcUaCommunicationConfig attribute*), 21  
 host (*TcpCommunicationConfig attribute*), 27  
 host (*TelnetCommunicationConfig attribute*), 29  
 host (*VisaCommunicationConfig attribute*), 32  
 hPascal (*PfeifferTPG.PressureUnits attribute*), 156  
 HT (*NewportConfigCommands attribute*), 142  
 hvl\_ccb  
     module, 191  
 hvl\_ccb.comm  
     module, 33  
 hvl\_ccb.comm.base  
     module, 11  
 hvl\_ccb.comm.labjack\_ljm  
     module, 15  
 hvl\_ccb.comm.modbus\_tcp  
     module, 18  
 hvl\_ccb.comm.opc  
     module, 20  
 hvl\_ccb.comm.serial  
     module, 23  
 hvl\_ccb.comm.tcp  
     module, 26  
 hvl\_ccb.comm.telnet  
     module, 28  
 hvl\_ccb.comm.visa  
     module, 30  
 hvl\_ccb.configuration  
     module, 187  
 hvl\_ccb.dev  
     module, 182  
 hvl\_ccb.dev.base  
     module, 92  
 hvl\_ccb.dev.crylas  
     module, 94  
 hvl\_ccb.dev.cube  
     module, 53  
 hvl\_ccb.dev.cube.base  
     module, 33  
 hvl\_ccb.dev.cube.constants  
     module, 38  
 hvl\_ccb.dev.cube.picube  
     module, 50  
 hvl\_ccb.dev.ea\_psi9000  
     module, 104  
 hvl\_ccb.dev.fug  
     module, 108  
 hvl\_ccb.dev.heinzinger  
     module, 119  
 hvl\_ccb.dev.highland\_t560  
     module, 58  
 hvl\_ccb.dev.highland\_t560.base  
     module, 53  
 hvl\_ccb.dev.highland\_t560.channel  
     module, 55  
 hvl\_ccb.dev.highland\_t560.device  
     module, 56  
 hvl\_ccb.dev.labjack  
     module, 125  
 hvl\_ccb.dev.lauda  
     module, 132  
 hvl\_ccb.dev.mbw973  
     module, 138  
 hvl\_ccb.dev.newport  
     module, 142  
 hvl\_ccb.dev.pfeiffer\_tpg  
     module, 155  
 hvl\_ccb.dev.rs\_rto1024  
     module, 161  
 hvl\_ccb.dev.se\_ils2t  
     module, 169  
 hvl\_ccb.dev.sst\_luminox  
     module, 175  
 hvl\_ccb.dev.supercube2015  
     module, 73  
 hvl\_ccb.dev.supercube2015.base  
     module, 58  
 hvl\_ccb.dev.supercube2015.constants  
     module, 64  
 hvl\_ccb.dev.supercube2015.typ\_a  
     module, 71  
 hvl\_ccb.dev.technix  
     module, 79  
 hvl\_ccb.dev.technix.base  
     module, 74  
 hvl\_ccb.dev.technix.device  
     module, 77  
 hvl\_ccb.dev.tiepie  
     module, 91  
 hvl\_ccb.dev.tiepie.base  
     module, 80  
 hvl\_ccb.dev.tiepie.channel  
     module, 82  
 hvl\_ccb.dev.tiepie.device  
     module, 84  
 hvl\_ccb.dev.tiepie.generator  
     module, 85  
 hvl\_ccb.dev.tiepie.i2c  
     module, 87  
 hvl\_ccb.dev.tiepie.oscilloscope  
     module, 88  
 hvl\_ccb.dev.tiepie.utils  
     module, 90  
 hvl\_ccb.dev.utils

- module, 179
  - `hvl_ccb.dev.visa`
    - module, 180
  - `hvl_ccb.exception`
    - module, 189
  - `hvl_ccb.experiment_manager`
    - module, 189
  - `hvl_ccb.utils`
    - module, 187
  - `hvl_ccb.utils.conversion_sensor`
    - module, 182
  - `hvl_ccb.utils.conversion_unit`
    - module, 183
  - `hvl_ccb.utils.enum`
    - module, 185
  - `hvl_ccb.utils.typing`
    - module, 186
  - `hvl_ccb.utils.validation`
    - module, 186
  - `hysteresis_compensation` (*NewportSMC100PPConfig attribute*), 150
- |
- I2C (*TiePieDeviceType attribute*), 81
  - ID (*FuGProbusIVCommands attribute*), 114
  - `Identification_error` (*PfeifferTPG.SensorStatus attribute*), 156
  - `identifier` (*LJMCommunicationConfig attribute*), 17
  - `identify_device()` (*FuG method*), 109
  - `identify_device()` (*HeinzingerPNC method*), 123
  - `identify_sensors()` (*PfeifferTPG method*), 157
  - IKR (*PfeifferTPG.SensorTypes attribute*), 156
  - IKR11 (*PfeifferTPG.SensorTypes attribute*), 156
  - IKR9 (*PfeifferTPG.SensorTypes attribute*), 156
  - ILS2T (*class in hvl\_ccb.dev.se\_ils2t*), 169
  - ILS2T.ActionsPtp (*class in hvl\_ccb.dev.se\_ils2t*), 169
  - ILS2T.Mode (*class in hvl\_ccb.dev.se\_ils2t*), 169
  - ILS2T.Ref16Jog (*class in hvl\_ccb.dev.se\_ils2t*), 169
  - ILS2T.State (*class in hvl\_ccb.dev.se\_ils2t*), 170
  - ILS2TConfig (*class in hvl\_ccb.dev.se\_ils2t*), 172
  - ILS2TException, 173
  - ILS2TModbusTcpCommunication (*class in hvl\_ccb.dev.se\_ils2t*), 173
  - ILS2TModbusTcpCommunicationConfig (*class in hvl\_ccb.dev.se\_ils2t*), 173
  - ILS2TRegAddr (*class in hvl\_ccb.dev.se\_ils2t*), 174
  - ILS2TRegDatatype (*class in hvl\_ccb.dev.se\_ils2t*), 174
  - IMMEDIATELY (*FuGRampModes attribute*), 117
  - IMPULSE\_140KV (*PowerSetup attribute*), 49
  - IMR (*PfeifferTPG.SensorTypes attribute*), 156
  - `in_1_1` (*GeneralSupport attribute*), 67
  - `in_1_2` (*GeneralSupport attribute*), 67
  - `in_2_1` (*GeneralSupport attribute*), 67
  - `in_2_2` (*GeneralSupport attribute*), 67
  - `in_3_1` (*GeneralSupport attribute*), 67
  - `in_3_2` (*GeneralSupport attribute*), 67
  - `in_4_1` (*GeneralSupport attribute*), 67
  - `in_4_2` (*GeneralSupport attribute*), 67
  - `in_5_1` (*GeneralSupport attribute*), 67
  - `in_5_2` (*GeneralSupport attribute*), 67
  - `in_6_1` (*GeneralSupport attribute*), 67
  - `in_6_2` (*GeneralSupport attribute*), 68
  - INACTIVE (*CryLasLaser.AnswersStatus attribute*), 97
  - INACTIVE (*DoorStatus attribute*), 47
  - inactive (*DoorStatus attribute*), 65
  - INACTIVE (*EarthingStickStatus attribute*), 48
  - inactive (*EarthingStickStatus attribute*), 67
  - inhibit (*Technix property*), 77
  - `init_attenuation` (*CryLasAttenuatorConfig attribute*), 95
  - `init_monitored_nodes()` (*OpcUaCommunication method*), 20
  - `init_shutter_status` (*CryLasLaserConfig attribute*), 101
  - `initialize()` (*NewportSMC100PP method*), 146
  - INITIALIZED (*ExperimentStatus attribute*), 190
  - INITIALIZING (*ExperimentStatus attribute*), 190
  - INITIALIZING (*SafetyStatus attribute*), 49
  - Initializing (*SafetyStatus attribute*), 70
  - INPUT (*FuGProbusVRegisterGroups attribute*), 116
  - INPUT (*GateMode attribute*), 53
  - `input()` (*GeneralSupport class method*), 68
  - `input_1` (*MeasurementsDividerRatio attribute*), 68
  - `input_1` (*MeasurementsScaledInput attribute*), 69
  - `input_2` (*MeasurementsScaledInput attribute*), 69
  - `input_3` (*MeasurementsScaledInput attribute*), 69
  - `input_4` (*MeasurementsScaledInput attribute*), 69
  - `input_range` (*TiePieOscilloscopeChannelConfig property*), 83
  - INSTALL (*AutoInstallMode attribute*), 53
  - INT32 (*ILS2TRegDatatype attribute*), 175
  - INT\_SYNTHESIZER (*TriggerMode attribute*), 55
  - `interface_type` (*PSI9000VisaCommunicationConfig attribute*), 107
  - `interface_type` (*RTO1024VisaCommunicationConfig attribute*), 168
  - `interface_type` (*VisaCommunicationConfig attribute*), 32
  - internal (*LabJack.CjcType attribute*), 126
  - INTERNAL (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 135
  - Internal (*PowerSetup attribute*), 70
  - InvalidSupercubeStatusError, 58
  - IO\_SCANNING (*ILS2TRegAddr attribute*), 174
  - IoScanningModeValueError, 175
  - `is_configdataclass` (*AsyncCommunicationProtocolConfig attribute*), 13

- is\_configdataclass (*BaseCubeConfiguration* attribute), 36  
 is\_configdataclass (*CryLasAttenuatorConfig* attribute), 95  
 is\_configdataclass (*CryLasLaserConfig* attribute), 101  
 is\_configdataclass (*EmptyConfig* attribute), 94, 188  
 is\_configdataclass (*FuGConfig* attribute), 110  
 is\_configdataclass (*HeinzingerConfig* attribute), 120  
 is\_configdataclass (*ILS2TConfig* attribute), 173  
 is\_configdataclass (*LaudaProRp245eConfig* attribute), 136  
 is\_configdataclass (*LJMCommunicationConfig* attribute), 17  
 is\_configdataclass (*LuminoxConfig* attribute), 177  
 is\_configdataclass (*MBW973Config* attribute), 140  
 is\_configdataclass (*ModbusTcpCommunicationConfig* attribute), 19  
 is\_configdataclass (*NewportSMC100PPConfig* attribute), 150  
 is\_configdataclass (*OpcUaCommunicationConfig* attribute), 21  
 is\_configdataclass (*PfeifferTPGConfig* attribute), 158  
 is\_configdataclass (*SupercubeConfiguration* attribute), 62  
 is\_configdataclass (*T560Config* attribute), 57  
 is\_configdataclass (*TcpCommunicationConfig* attribute), 27  
 is\_configdataclass (*TechnixConfig* attribute), 78  
 is\_configdataclass (*TiePieDeviceConfig* attribute), 80  
 is\_configdataclass (*VisaCommunicationConfig* attribute), 32  
 is\_data\_ready\_polling\_interval\_sec (*TiePieDeviceConfig* attribute), 80  
 is\_done() (*MBW973* method), 138  
 is\_error() (*ExperimentManager* method), 190  
 is\_finished() (*ExperimentManager* method), 190  
 is\_generic\_type\_hint() (in module *hvl\_ccb.utils*.typing), 186  
 is\_in\_range() (*ILS2TRegDatatype* method), 175  
 is\_inactive (*CryLasLaser.LaserStatus* property), 98  
 is\_measurement\_data\_ready() (*TiePieOscilloscope* method), 89  
 is\_open (*Client* property), 20  
 is\_open (*LJMCommunication* property), 15  
 is\_open (*OpcUaCommunication* property), 20  
 is\_open (*SerialCommunication* property), 23  
 is\_open (*TelnetCommunication* property), 28  
 is\_polling() (*Poller* method), 179  
 is\_ready (*CryLasLaser.LaserStatus* property), 98  
 is\_running() (*ExperimentManager* method), 190  
 is\_started (*Technix* property), 77  
 is\_triggered() (*TiePieOscilloscope* method), 89  
 is\_valid\_scale\_range\_reversed\_str() (*PfeifferTPGConfig.Model* method), 158
- ## J
- J (*LabJack.ThermocoupleType* attribute), 128  
 jerk\_time (*NewportSMC100PPConfig* attribute), 150  
 JOG (*ILS2T.Mode* attribute), 169  
 jog\_run() (*ILS2T* method), 171  
 jog\_stop() (*ILS2T* method), 171  
 JOGGING (*NewportStates* attribute), 155  
 JOGGING\_FROM\_DISABLE (*NewportSMC100PP.StateMessages* attribute), 143  
 JOGGING\_FROM\_READY (*NewportSMC100PP.StateMessages* attribute), 143  
 JOGN\_FAST (*ILS2TRegAddr* attribute), 174  
 JOGN\_SLOW (*ILS2TRegAddr* attribute), 174  
 JR (*NewportConfigCommands* attribute), 142
- ## K
- K (*LabJack.TemperatureUnit* attribute), 127  
 K (*LabJack.ThermocoupleType* attribute), 128  
 K (*Temperature* attribute), 184  
 KELVIN (*Temperature* attribute), 184  
 keys() (*AsyncCommunicationProtocolConfig* class method), 13  
 keys() (*BaseCubeConfiguration* class method), 36  
 keys() (*BaseCubeOpcUaCommunicationConfig* class method), 37  
 keys() (*CryLasAttenuatorConfig* class method), 95  
 keys() (*CryLasAttenuatorSerialCommunicationConfig* class method), 97  
 keys() (*CryLasLaserConfig* class method), 101  
 keys() (*CryLasLaserSerialCommunicationConfig* class method), 103  
 keys() (*EmptyConfig* class method), 94, 188  
 keys() (*FuGConfig* class method), 110  
 keys() (*FuGSerialCommunicationConfig* class method), 118  
 keys() (*HeinzingerConfig* class method), 120  
 keys() (*HeinzingerSerialCommunicationConfig* class method), 124  
 keys() (*ILS2TConfig* class method), 173  
 keys() (*ILS2TModbusTcpCommunicationConfig* class method), 174  
 keys() (*LaudaProRp245eConfig* class method), 136  
 keys() (*LaudaProRp245eTcpCommunicationConfig* class method), 137  
 keys() (*LJMCommunicationConfig* class method), 17  
 keys() (*LuminoxConfig* class method), 177  
 keys() (*LuminoxSerialCommunicationConfig* class method), 179

- keys() (*MBW973Config* class method), 140
- keys() (*MBW973SerialCommunicationConfig* class method), 141
- keys() (*ModbusTcpCommunicationConfig* class method), 19
- keys() (*NewportSMC100PPConfig* class method), 150
- keys() (*NewportSMC100PPSerialCommunicationConfig* class method), 154
- keys() (*OpCuaCommunicationConfig* class method), 21
- keys() (*PfeifferTPGConfig* class method), 158
- keys() (*PfeifferTPGSerialCommunicationConfig* class method), 160
- keys() (*PICubeConfiguration* class method), 51
- keys() (*PICubeOpCuaCommunicationConfig* class method), 52
- keys() (*PSI9000Config* class method), 106
- keys() (*PSI9000VisaCommunicationConfig* class method), 108
- keys() (*RTO1024Config* class method), 167
- keys() (*RTO1024VisaCommunicationConfig* class method), 168
- keys() (*SerialCommunicationConfig* class method), 25
- keys() (*SupercubeAOpCuaConfiguration* class method), 73
- keys() (*SupercubeConfiguration* class method), 62
- keys() (*SupercubeOpCuaCommunicationConfig* class method), 63
- keys() (*T560CommunicationConfig* class method), 54
- keys() (*T560Config* class method), 57
- keys() (*TcpCommunicationConfig* class method), 27
- keys() (*TechnixConfig* class method), 78
- keys() (*TechnixSerialCommunicationConfig* class method), 75
- keys() (*TechnixTelnetCommunicationConfig* class method), 76
- keys() (*TelnetCommunicationConfig* class method), 29
- keys() (*TiePieDeviceConfig* class method), 80
- keys() (*VisaCommunicationConfig* class method), 32
- keys() (*VisaDeviceConfig* class method), 182
- kV (*HeinzingerPNC.UnitVoltage* attribute), 123
- L**
- LabJack (class in *hvl\_ccb.dev.labjack*), 126
- LabJack.AInRange (class in *hvl\_ccb.dev.labjack*), 126
- LabJack.BitLimit (class in *hvl\_ccb.dev.labjack*), 126
- LabJack.CalMicroAmpere (class in *hvl\_ccb.dev.labjack*), 126
- LabJack.CjcType (class in *hvl\_ccb.dev.labjack*), 126
- LabJack.ClockFrequency (class in *hvl\_ccb.dev.labjack*), 126
- LabJack.DeviceType (class in *hvl\_ccb.dev.labjack*), 127
- LabJack.DIOStatus (class in *hvl\_ccb.dev.labjack*), 127
- LabJack.TemperatureUnit (class in *hvl\_ccb.dev.labjack*), 127
- LabJack.ThermocoupleType (class in *hvl\_ccb.dev.labjack*), 127
- LabJackError, 131
- LabJackIdentifierDIOError, 131
- laser\_off() (*CryLasLaser* method), 98
- laser\_on() (*CryLasLaser* method), 99
- LaudaProRp245e (class in *hvl\_ccb.dev.lauda*), 132
- LaudaProRp245eCommand (class in *hvl\_ccb.dev.lauda*), 134
- LaudaProRp245eCommandError, 135
- LaudaProRp245eConfig (class in *hvl\_ccb.dev.lauda*), 135
- LaudaProRp245eConfig.ExtControlModeEnum (class in *hvl\_ccb.dev.lauda*), 135
- LaudaProRp245eConfig.OperationModeEnum (class in *hvl\_ccb.dev.lauda*), 136
- LaudaProRp245eTcpCommunication (class in *hvl\_ccb.dev.lauda*), 137
- LaudaProRp245eTcpCommunicationConfig (class in *hvl\_ccb.dev.lauda*), 137
- LEM4000S (class in *hvl\_ccb.utils.conversion\_sensor*), 182
- LF (*FuGTerminators* attribute), 119
- LFCR (*FuGTerminators* attribute), 119
- LINE\_1 (*MessageBoard* attribute), 48
- LINE\_10 (*MessageBoard* attribute), 48
- LINE\_11 (*MessageBoard* attribute), 48
- LINE\_12 (*MessageBoard* attribute), 48
- LINE\_13 (*MessageBoard* attribute), 48
- LINE\_14 (*MessageBoard* attribute), 48
- LINE\_15 (*MessageBoard* attribute), 48
- LINE\_2 (*MessageBoard* attribute), 48
- LINE\_3 (*MessageBoard* attribute), 48
- LINE\_4 (*MessageBoard* attribute), 48
- LINE\_5 (*MessageBoard* attribute), 48
- LINE\_6 (*MessageBoard* attribute), 48
- LINE\_7 (*MessageBoard* attribute), 48
- LINE\_8 (*MessageBoard* attribute), 48
- LINE\_9 (*MessageBoard* attribute), 48
- list\_devices() (*TiePieOscilloscope* static method), 89
- list\_directory() (*RTO1024* method), 163
- LJMCommunication (class in *hvl\_ccb.comm.labjack\_ljm*), 15
- LJMCommunicationConfig (class in *hvl\_ccb.comm.labjack\_ljm*), 16
- LJMCommunicationConfig.ConnectionType (class in *hvl\_ccb.comm.labjack\_ljm*), 16
- LJMCommunicationConfig.DeviceType (class in *hvl\_ccb.comm.labjack\_ljm*), 16
- LJMCommunicationError, 17
- lm34 (*LabJack.CjcType* attribute), 126
- LMT70A (class in *hvl\_ccb.utils.conversion\_sensor*), 183
- load\_configuration() (*RTO1024* method), 164

- load\_device\_configuration() (*T560 method*), 57
  - local\_display() (*RTO1024 method*), 164
  - LOCKED (*DoorStatus attribute*), 47
  - locked (*DoorStatus attribute*), 65
  - LOW (*LabJack.DIOStatus attribute*), 127
  - LOWER\_TEMP (*LaudaProRp245eCommand attribute*), 134
  - lower\_temp (*LaudaProRp245eConfig attribute*), 136
  - Luminox (*class in hvl\_ccb.dev.sst\_luminox*), 175
  - LuminoxConfig (*class in hvl\_ccb.dev.sst\_luminox*), 176
  - LuminoxException, 177
  - LuminoxMeasurementType (*class in hvl\_ccb.dev.sst\_luminox*), 177
  - LuminoxMeasurementTypeDict (*in module hvl\_ccb.dev.sst\_luminox*), 178
  - LuminoxMeasurementTypeError, 178
  - LuminoxMeasurementTypeValue (*in module hvl\_ccb.dev.sst\_luminox*), 178
  - LuminoxOutputMode (*class in hvl\_ccb.dev.sst\_luminox*), 178
  - LuminoxOutputModeError, 178
  - LuminoxSerialCommunication (*class in hvl\_ccb.dev.sst\_luminox*), 178
  - LuminoxSerialCommunicationConfig (*class in hvl\_ccb.dev.sst\_luminox*), 178
  - LUT (*LMT70A attribute*), 183
- ## M
- mA (*HeinzingerPNC.UnitCurrent attribute*), 122
  - MANUAL (*EarthingStickOperatingStatus attribute*), 47
  - manual() (*EarthingStick class method*), 65
  - manual\_1 (*EarthingStick attribute*), 65
  - manual\_2 (*EarthingStick attribute*), 65
  - manual\_3 (*EarthingStick attribute*), 65
  - manual\_4 (*EarthingStick attribute*), 65
  - manual\_5 (*EarthingStick attribute*), 65
  - manual\_6 (*EarthingStick attribute*), 66
  - MARK (*SerialCommunicationParity attribute*), 25
  - max\_current (*FuG property*), 109
  - max\_current (*HeinzingerPNC property*), 123
  - max\_current (*Technix property*), 77
  - max\_current (*TechnixConfig attribute*), 79
  - max\_current\_hardware (*FuG property*), 109
  - max\_current\_hardware (*HeinzingerPNC property*), 123
  - max\_pr\_number (*LaudaProRp245eConfig attribute*), 136
  - max\_pump\_level (*LaudaProRp245eConfig attribute*), 136
  - max\_timeout\_retry\_nr (*OpcUaCommunicationConfig attribute*), 22
  - max\_voltage (*FuG property*), 109
  - max\_voltage (*HeinzingerPNC property*), 123
  - max\_voltage (*Technix property*), 77
  - max\_voltage (*TechnixConfig attribute*), 79
  - max\_voltage\_hardware (*FuG property*), 109
  - max\_voltage\_hardware (*HeinzingerPNC property*), 123
  - MAXIMUM (*LabJack.ClockFrequency attribute*), 127
  - mbar (*PfeifferTPG.PressureUnits attribute*), 156
  - MBW973 (*class in hvl\_ccb.dev.mbw973*), 138
  - MBW973Config (*class in hvl\_ccb.dev.mbw973*), 139
  - MBW973ControlRunningException, 140
  - MBW973Error, 140
  - MBW973PumpRunningException, 140
  - MBW973SerialCommunication (*class in hvl\_ccb.dev.mbw973*), 140
  - MBW973SerialCommunicationConfig (*class in hvl\_ccb.dev.mbw973*), 140
  - measure() (*PfeifferTPG method*), 157
  - measure\_all() (*PfeifferTPG method*), 157
  - measure\_current() (*HeinzingerDI method*), 121
  - measure\_voltage() (*HeinzingerDI method*), 121
  - measure\_voltage\_current() (*PSI9000 method*), 105
  - MeasurementsDividerRatio (*class in hvl\_ccb.dev.supercube2015.constants*), 68
  - MeasurementsScaledInput (*class in hvl\_ccb.dev.supercube2015.constants*), 68
  - MessageBoard (*class in hvl\_ccb.dev.cube.constants*), 48
  - micro\_step\_per\_full\_step\_factor (*NewportSMC100PPConfig attribute*), 150
  - Micron (*PfeifferTPG.PressureUnits attribute*), 155
  - MILLIMETER\_MERCURY (*Pressure attribute*), 184
  - MINIMUM (*LabJack.ClockFrequency attribute*), 127
  - MMHG (*Pressure attribute*), 184
  - ModbusTcpCommunication (*class in hvl\_ccb.comm.modbus\_tcp*), 18
  - ModbusTcpCommunicationConfig (*class in hvl\_ccb.comm.modbus\_tcp*), 19
  - ModbusTcpConnectionFailedException, 19
  - model (*PfeifferTPGConfig attribute*), 159
  - module
    - hvl\_ccb, 191
    - hvl\_ccb.comm, 33
    - hvl\_ccb.comm.base, 11
    - hvl\_ccb.comm.labjack\_ljm, 15
    - hvl\_ccb.comm.modbus\_tcp, 18
    - hvl\_ccb.comm.opc, 20
    - hvl\_ccb.comm.serial, 23
    - hvl\_ccb.comm.tcp, 26
    - hvl\_ccb.comm.telnet, 28
    - hvl\_ccb.comm.visa, 30
    - hvl\_ccb.configuration, 187
    - hvl\_ccb.dev, 182
    - hvl\_ccb.dev.base, 92
    - hvl\_ccb.dev.crylas, 94
    - hvl\_ccb.dev.cube, 53
    - hvl\_ccb.dev.cube.base, 33
    - hvl\_ccb.dev.cube.constants, 38
    - hvl\_ccb.dev.cube.picube, 50

- hvl\_ccb.dev.ea\_psi9000, 104
  - hvl\_ccb.dev.fug, 108
  - hvl\_ccb.dev.heinzinger, 119
  - hvl\_ccb.dev.highland\_t560, 58
  - hvl\_ccb.dev.highland\_t560.base, 53
  - hvl\_ccb.dev.highland\_t560.channel, 55
  - hvl\_ccb.dev.highland\_t560.device, 56
  - hvl\_ccb.dev.labjack, 125
  - hvl\_ccb.dev.lauda, 132
  - hvl\_ccb.dev.mbw973, 138
  - hvl\_ccb.dev.newport, 142
  - hvl\_ccb.dev.pfeiffer\_tpg, 155
  - hvl\_ccb.dev.rs\_rto1024, 161
  - hvl\_ccb.dev.se\_ils2t, 169
  - hvl\_ccb.dev.sst\_luminox, 175
  - hvl\_ccb.dev.supercube2015, 73
  - hvl\_ccb.dev.supercube2015.base, 58
  - hvl\_ccb.dev.supercube2015.constants, 64
  - hvl\_ccb.dev.supercube2015.typ\_a, 71
  - hvl\_ccb.dev.technix, 79
  - hvl\_ccb.dev.technix.base, 74
  - hvl\_ccb.dev.technix.device, 77
  - hvl\_ccb.dev.tiepie, 91
  - hvl\_ccb.dev.tiepie.base, 80
  - hvl\_ccb.dev.tiepie.channel, 82
  - hvl\_ccb.dev.tiepie.device, 84
  - hvl\_ccb.dev.tiepie.generator, 85
  - hvl\_ccb.dev.tiepie.i2c, 87
  - hvl\_ccb.dev.tiepie.oscilloscope, 88
  - hvl\_ccb.dev.tiepie.utils, 90
  - hvl\_ccb.dev.utils, 179
  - hvl\_ccb.dev.visa, 180
  - hvl\_ccb.exception, 189
  - hvl\_ccb.experiment\_manager, 189
  - hvl\_ccb.utils, 187
  - hvl\_ccb.utils.conversion\_sensor, 182
  - hvl\_ccb.utils.conversion\_unit, 183
  - hvl\_ccb.utils.enum, 185
  - hvl\_ccb.utils.typing, 186
  - hvl\_ccb.utils.validation, 186
  - MONITOR\_I (*FuGProbusVRegisterGroups* attribute), 116
  - MONITOR\_V (*FuGProbusVRegisterGroups* attribute), 116
  - most\_recent\_error (*FuGProbusVConfigRegisters* property), 114
  - motion\_distance\_per\_full\_step (*NewportSMC100PPConfig* attribute), 150
  - motor\_config (*NewportSMC100PPConfig* property), 150
  - move\_to\_absolute\_position() (*NewportSMC100PP* method), 147
  - move\_to\_relative\_position() (*NewportSMC100PP* method), 147
  - move\_wait\_sec (*NewportSMC100PPConfig* attribute), 150
  - MOVING (*NewportSMC100PP.StateMessages* attribute), 143
  - MOVING (*NewportStates* attribute), 155
  - MS\_NOMINAL\_CURRENT (*PSI9000* attribute), 104
  - MS\_NOMINAL\_VOLTAGE (*PSI9000* attribute), 104
  - MULTI\_COMMANDS\_MAX (*VisaCommunication* attribute), 30
  - MULTI\_COMMANDS\_SEPARATOR (*VisaCommunication* attribute), 30
- ## N
- n\_channels (*TiePieOscilloscope* property), 89
  - n\_max\_try\_get\_device (*TiePieDeviceConfig* attribute), 80
  - NameEnum (*class in hvl\_ccb.utils.enum*), 185
  - NAMES (*SerialCommunicationParity* attribute), 25
  - names() (*RTO1024.TriggerModes* class method), 162
  - namespace\_index (*BaseCubeConfiguration* attribute), 36
  - namespace\_index (*SupercubeConfiguration* attribute), 62
  - NATIVEONLY (*TiePieOscilloscopeAutoResolutionModes* attribute), 89
  - NED\_END\_OF\_TURN (*NewportSMC100PP.MotorErrors* attribute), 143
  - NEG (*ILS2T.Ref16Jog* attribute), 169
  - NEG\_FAST (*ILS2T.Ref16Jog* attribute), 169
  - NEGATIVE (*FuGPolarities* attribute), 113
  - NEGATIVE (*Polarity* attribute), 48
  - negative\_software\_limit (*NewportSMC100PPConfig* attribute), 150
  - NewportConfigCommands (*class in hvl\_ccb.dev.newport*), 142
  - NewportControllerError, 142
  - NewportError, 142
  - NewportMotorError, 142
  - NewportMotorPowerSupplyWasCutError, 143
  - NewportSerialCommunicationError, 154
  - NewportSMC100PP (*class in hvl\_ccb.dev.newport*), 143
  - NewportSMC100PP.MotorErrors (*class in hvl\_ccb.dev.newport*), 143
  - NewportSMC100PP.StateMessages (*class in hvl\_ccb.dev.newport*), 143
  - NewportSMC100PPConfig (*class in hvl\_ccb.dev.newport*), 149
  - NewportSMC100PPConfig.EspStageConfig (*class in hvl\_ccb.dev.newport*), 149
  - NewportSMC100PPConfig.HomeSearch (*class in hvl\_ccb.dev.newport*), 149
  - NewportSMC100PPSerialCommunication (*class in hvl\_ccb.dev.newport*), 151
  - NewportSMC100PPSerialCommunication.ControllerErrors (*class in hvl\_ccb.dev.newport*), 151

NewportSMC100PPSerialCommunicationConfig (class in *hvl\_ccb.dev.newport*), 153

NewportStates (class in *hvl\_ccb.dev.newport*), 154

NewportUncertainPositionError, 155

NO (*FuGDigitalVal* attribute), 111

NO\_ERROR (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 151

NO\_REF (*NewportStates* attribute), 155

NO\_REF\_ESP\_STAGE\_ERROR (New-attribute), 143

NO\_REF\_FROM\_CONFIG (New-attribute), 143

NO\_REF\_FROM\_DISABLED (New-attribute), 143

NO\_REF\_FROM\_HOMING (New-attribute), 144

NO\_REF\_FROM\_JOGGING (New-attribute), 144

NO\_REF\_FROM\_MOVING (New-attribute), 144

NO\_REF\_FROM\_READY (New-attribute), 144

NO\_REF\_FROM\_RESET (New-attribute), 144

No\_sensor (*PfeifferTPG.SensorStatus* attribute), 156

NO\_SOURCE (*PowerSetup* attribute), 49

NOISE (*TiePieGeneratorSignalType* attribute), 86

noise\_level\_measurement\_channel\_1 (*BaseCube-Configuration* attribute), 36

noise\_level\_measurement\_channel\_2 (*BaseCube-Configuration* attribute), 36

noise\_level\_measurement\_channel\_3 (*BaseCube-Configuration* attribute), 36

noise\_level\_measurement\_channel\_4 (*BaseCube-Configuration* attribute), 36

NONE (*ILS2T.Ref16Jog* attribute), 169

NONE (*LabJack.ThermocoupleType* attribute), 128

None (*PfeifferTPG.SensorTypes* attribute), 156

NONE (*SerialCommunicationParity* attribute), 25

NORMAL (*RTO1024.TriggerModes* attribute), 162

noSen (*PfeifferTPG.SensorTypes* attribute), 156

noSENSOR (*PfeifferTPG.SensorTypes* attribute), 156

not\_defined (*AlarmText* attribute), 65

nr\_trials\_activate (*LuminoxConfig* attribute), 177

NullCommunicationProtocol (class in *hvl\_ccb.comm.base*), 14

Number (in module *hvl\_ccb.utils.typing*), 186

number\_of\_decimals (*HeinzingerConfig* attribute), 120

number\_of\_sensors (*PfeifferTPG* property), 157

**O**

OFF (*SerialCommunicationParity* attribute), 25

OFF (*AutoInstallMode* attribute), 53

OFF (*FuGDigitalVal* attribute), 111

OFF (*GateMode* attribute), 53

OFF (*HeinzingerDI.OutputStatus* attribute), 121

OFF (*TriggerMode* attribute), 55

offset (*TiePieGeneratorConfig* property), 86

OH (*NewportConfigCommands* attribute), 142

Ok (*PfeifferTPG.SensorStatus* attribute), 156

on (*FuG* property), 109

ON (*FuGDigitalVal* attribute), 111

on (*FuGProbusVDIRegisters* property), 115

ON (*HeinzingerDI.OutputStatus* attribute), 121

ON (*ILS2T.State* attribute), 170

ONE (*HeinzingerConfig.RecordingsEnum* attribute), 120

ONE (*LabJack.AInRange* attribute), 126

ONE (*SerialCommunicationStopbits* attribute), 26

ONE\_HUNDREDTH (*LabJack.AInRange* attribute), 126

ONE\_POINT\_FIVE (*SerialCommunicationStopbits* attribute), 26

ONE\_TENTH (*LabJack.AInRange* attribute), 126

ONLYUPWARDSOFFTOZERO (*FuGRampModes* attribute), 117

OPC\_MAX\_YEAR (*BaseCube* attribute), 33

OPC\_MIN\_YEAR (*BaseCube* attribute), 33

OpcUaCommunication (class in *hvl\_ccb.comm.opc*), 20

OpcUaCommunicationConfig (class in *hvl\_ccb.comm.opc*), 21

OpcUaCommunicationIOError, 22

OpcUaCommunicationTimeoutError, 22

OpcUaSubHandler (class in *hvl\_ccb.comm.opc*), 22

OPEN (*DoorStatus* attribute), 47

open (*DoorStatus* attribute), 65

OPEN (*EarthingStickOperation* attribute), 47

OPEN (*EarthingStickStatus* attribute), 48

open (*EarthingStickStatus* attribute), 67

open() (*CommunicationProtocol* method), 14

open() (*LaudaProRp245eTcpCommunication* method), 137

open() (*LJMCommunication* method), 15

open() (*ModbusTcpCommunication* method), 18

open() (*NullCommunicationProtocol* method), 14

open() (*OpcUaCommunication* method), 20

open() (*SerialCommunication* method), 23

open() (*Tcp* method), 26

open() (*TelnetCommunication* method), 28

open() (*VisaCommunication* method), 30

open\_interlock (*Technix* property), 78

open\_shutter() (*CryLasLaser* method), 99

- `open_timeout` (*VisaCommunicationConfig* attribute), 32
- `OPENED` (*CryLasLaser.AnswersShutter* attribute), 97
- `OPENED` (*CryLasLaserShutterStatus* attribute), 103
- `operate` (*BaseCube* property), 35
- `operate` (*PICube* property), 50
- `operate()` (*Supercube2015Base* method), 60
- `OPERATION_MODE` (*LaudaProRp245eCommand* attribute), 134
- `operation_mode` (*LaudaProRp245eConfig* attribute), 136
- `optional_defaults()` (*AsyncCommunicationProtocolConfig* class method), 13
- `optional_defaults()` (*BaseCubeConfiguration* class method), 36
- `optional_defaults()` (*BaseCubeOpcUaCommunicationConfig* class method), 37
- `optional_defaults()` (*CryLasAttenuatorConfig* class method), 96
- `optional_defaults()` (*CryLasAttenuatorSerialCommunicationConfig* class method), 97
- `optional_defaults()` (*CryLasLaserConfig* class method), 101
- `optional_defaults()` (*CryLasLaserSerialCommunicationConfig* class method), 103
- `optional_defaults()` (*EmptyConfig* class method), 94, 188
- `optional_defaults()` (*FuGConfig* class method), 110
- `optional_defaults()` (*FuGSerialCommunicationConfig* class method), 118
- `optional_defaults()` (*HeinzingerConfig* class method), 120
- `optional_defaults()` (*HeinzingerSerialCommunicationConfig* class method), 124
- `optional_defaults()` (*ILS2TConfig* class method), 173
- `optional_defaults()` (*ILS2TModbusTcpCommunicationConfig* class method), 174
- `optional_defaults()` (*LaudaProRp245eConfig* class method), 136
- `optional_defaults()` (*LaudaProRp245eTcpCommunicationConfig* class method), 138
- `optional_defaults()` (*LJMCommunicationConfig* class method), 17
- `optional_defaults()` (*LuminosConfig* class method), 177
- `optional_defaults()` (*LuminosSerialCommunicationConfig* class method), 179
- `optional_defaults()` (*MBW973Config* class method), 140
- `optional_defaults()` (*MBW973SerialCommunicationConfig* class method), 141
- `optional_defaults()` (*ModbusTcpCommunicationConfig* class method), 19
- `optional_defaults()` (*NewportSMC100PPConfig* class method), 150
- `optional_defaults()` (*NewportSMC100PPSerialCommunicationConfig* class method), 154
- `optional_defaults()` (*OpcUaCommunicationConfig* class method), 22
- `optional_defaults()` (*PfeifferTPGConfig* class method), 159
- `optional_defaults()` (*PfeifferTPGSerialCommunicationConfig* class method), 160
- `optional_defaults()` (*PICubeConfiguration* class method), 51
- `optional_defaults()` (*PICubeOpcUaCommunicationConfig* class method), 52
- `optional_defaults()` (*PSI9000Config* class method), 106
- `optional_defaults()` (*PSI9000VisaCommunicationConfig* class method), 108
- `optional_defaults()` (*RTO1024Config* class method), 167
- `optional_defaults()` (*RTO1024VisaCommunicationConfig* class method), 168
- `optional_defaults()` (*SerialCommunicationConfig* class method), 25
- `optional_defaults()` (*SupercubeAOpcUaConfiguration* class method), 73
- `optional_defaults()` (*SupercubeConfiguration* class method), 62
- `optional_defaults()` (*SupercubeOpcUaCommunicationConfig* class method), 63
- `optional_defaults()` (*T560CommunicationConfig* class method), 54
- `optional_defaults()` (*T560Config* class method), 57
- `optional_defaults()` (*TcpCommunicationConfig* class method), 27
- `optional_defaults()` (*TechnixConfig* class method), 79
- `optional_defaults()` (*TechnixSerialCommunicationConfig* class method), 75
- `optional_defaults()` (*TechnixTelnetCommunicationConfig* class method), 76
- `optional_defaults()` (*TelnetCommunicationConfig* class method), 29
- `optional_defaults()` (*TiePieDeviceConfig* class method), 81
- `optional_defaults()` (*VisaCommunicationConfig* class method), 32
- `optional_defaults()` (*VisaDeviceConfig* class method), 182

- OSCILLOSCOPE (*TiePieDeviceType* attribute), 81  
 OT (*NewportConfigCommands* attribute), 142  
 out (*FuGProbusVDORegisters* property), 115  
 out\_1\_1 (*GeneralSupport* attribute), 68  
 out\_1\_2 (*GeneralSupport* attribute), 68  
 out\_2\_1 (*GeneralSupport* attribute), 68  
 out\_2\_2 (*GeneralSupport* attribute), 68  
 out\_3\_1 (*GeneralSupport* attribute), 68  
 out\_3\_2 (*GeneralSupport* attribute), 68  
 out\_4\_1 (*GeneralSupport* attribute), 68  
 out\_4\_2 (*GeneralSupport* attribute), 68  
 out\_5\_1 (*GeneralSupport* attribute), 68  
 out\_5\_2 (*GeneralSupport* attribute), 68  
 out\_6\_1 (*GeneralSupport* attribute), 68  
 out\_6\_2 (*GeneralSupport* attribute), 68  
 OUTPUT (*FuGProbusIVCommands* attribute), 114  
 OUTPUT (*GateMode* attribute), 53  
 output (*Technix* property), 78  
 output() (*GeneralSupport* class method), 68  
 output\_off() (*FuGProbusIV* method), 113  
 output\_off() (*HeinzingerDI* method), 121  
 output\_on() (*HeinzingerDI* method), 122  
 OUTPUT\_POWER\_EXCEEDED (*NewportSMC100PP.MotorErrors* attribute), 143  
 output\_status (*HeinzingerDI* property), 122  
 OUTPUTONCMD (*FuGProbusVRegisterGroups* attribute), 116  
 OUTPUTX0 (*FuGProbusVRegisterGroups* attribute), 116  
 OUTPUTX1 (*FuGProbusVRegisterGroups* attribute), 116  
 OUTPUTX2 (*FuGProbusVRegisterGroups* attribute), 116  
 OUTPUTXCMD (*FuGProbusVRegisterGroups* attribute), 116  
 outX0 (*FuG* property), 109  
 outX1 (*FuG* property), 109  
 outX2 (*FuG* property), 109  
 outXCMD (*FuG* property), 109  
 Overrange (*PfeifferTPG.SensorStatus* attribute), 156
- P**
- PA (*Pressure* attribute), 184  
 PARAM\_MISSING\_OR\_INVALID (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 151  
 parity (*CryLasAttenuatorSerialCommunicationConfig* attribute), 97  
 parity (*CryLasLaserSerialCommunicationConfig* attribute), 103  
 parity (*FuGSerialCommunicationConfig* attribute), 118  
 parity (*HeinzingerSerialCommunicationConfig* attribute), 124  
 parity (*LuminoxSerialCommunicationConfig* attribute), 179  
 parity (*MBW973SerialCommunicationConfig* attribute), 141  
 parity (*NewportSMC100PPSerialCommunicationConfig* attribute), 154  
 parity (*PfeifferTPGSerialCommunicationConfig* attribute), 160  
 Parity (*SerialCommunicationConfig* attribute), 24  
 parity (*SerialCommunicationConfig* attribute), 25  
 parse\_read\_measurement\_value() (*LuminoxMeasurementType* method), 177  
 partial\_pressure\_o2 (*LuminoxMeasurementType* attribute), 177  
 Pascal (*PfeifferTPG.PressureUnits* attribute), 155  
 PASCAL (*Pressure* attribute), 184  
 pause() (*LaudaProRp245e* method), 132  
 pause\_ramp() (*LaudaProRp245e* method), 132  
 PBR (*PfeifferTPG.SensorTypes* attribute), 156  
 PEAK\_CURRENT\_LIMIT (*NewportSMC100PP.MotorErrors* attribute), 143  
 peak\_output\_current\_limit (*NewportSMC100PPConfig* attribute), 150  
 percent\_o2 (*LuminoxMeasurementType* attribute), 177  
 period (*T560* property), 57  
 PfeifferTPG (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 155  
 PfeifferTPG.PressureUnits (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 155  
 PfeifferTPG.SensorStatus (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 156  
 PfeifferTPG.SensorTypes (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 156  
 PfeifferTPGConfig (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 158  
 PfeifferTPGConfig.Model (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 158  
 PfeifferTPGError, 159  
 PfeifferTPGSerialCommunication (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 159  
 PfeifferTPGSerialCommunicationConfig (*class in hvl\_ccb.dev.pfeiffer\_tpg*), 159  
 PICube (*class in hvl\_ccb.dev.cube.picube*), 50  
 PICubeConfiguration (*class in hvl\_ccb.dev.cube.picube*), 51  
 PICubeOpcUaCommunication (*class in hvl\_ccb.dev.cube.picube*), 51  
 PICubeOpcUaCommunicationConfig (*class in hvl\_ccb.dev.cube.picube*), 52  
 PICubeTestParameterError, 48  
 PKR (*PfeifferTPG.SensorTypes* attribute), 156  
 Polarity (*class in hvl\_ccb.dev.cube.constants*), 48  
 Polarity (*class in hvl\_ccb.dev.highland\_t560.base*), 53  
 POLARITY (*FuGProbusIVCommands* attribute), 114  
 polarity (*PICube* property), 50  
 Poller (*class in hvl\_ccb.dev.utils*), 179

- polling (*LuminoxOutputMode* attribute), 178
- polling\_delay\_sec (*BaseCubeConfiguration* attribute), 37
- polling\_interval (*MBW973Config* attribute), 140
- polling\_interval\_sec (*BaseCubeConfiguration* attribute), 37
- polling\_interval\_sec (*TechnixConfig* attribute), 79
- polling\_period (*CryLasLaserConfig* attribute), 101
- polling\_timeout (*CryLasLaserConfig* attribute), 101
- port (*ModbusTcpCommunicationConfig* attribute), 19
- port (*OpcUaCommunicationConfig* attribute), 22
- port (*SerialCommunicationConfig* attribute), 25
- port (*SupercubeOpcUaCommunicationConfig* attribute), 63
- port (*T560CommunicationConfig* attribute), 54
- port (*TcpCommunicationConfig* attribute), 27
- port (*TechnixTelnetCommunicationConfig* attribute), 76
- port (*TelnetCommunicationConfig* attribute), 29
- port (*VisaCommunicationConfig* attribute), 32
- POS (*ILS2T.Ref16Jog* attribute), 170
- POS\_END\_OF\_TURN (*NewportSMC100PP.MotorErrors* attribute), 143
- POS\_FAST (*ILS2T.Ref16Jog* attribute), 170
- POSITION (*ILS2TRegAddr* attribute), 174
- POSITION\_OUT\_OF\_LIMIT (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 151
- POSITIVE (*FuGPolarities* attribute), 113
- POSITIVE (*Polarity* attribute), 48
- positive\_software\_limit (*NewportSMC100PPConfig* attribute), 150
- post\_force\_value() (*NewportSMC100PPConfig* method), 150
- post\_stop\_pause\_sec (*TechnixConfig* attribute), 79
- POUNDS\_PER\_SQUARE\_INCH (*Pressure* attribute), 184
- Power (*class in hvl\_ccb.dev.supercube2015.constants*), 69
- POWER\_INVERTER\_220V (*PowerSetup* attribute), 49
- power\_limit (*PSI9000Config* attribute), 107
- power\_setup (*PICube* property), 51
- PowerSetup (*class in hvl\_ccb.dev.cube.constants*), 48
- PowerSetup (*class in hvl\_ccb.dev.supercube2015.constants*), 69
- pre\_sample\_ratio (*TiePieOscilloscopeConfig* property), 90
- prepare\_ultra\_segmentation() (*RTO1024* method), 164
- preserve\_type() (*in module hvl\_ccb.utils.conversion\_unit*), 184
- Pressure (*class in hvl\_ccb.utils.conversion\_unit*), 183
- probe\_offset (*TiePieOscilloscopeChannelConfig* property), 83
- PSI (*Pressure* attribute), 184
- PSI9000 (*class in hvl\_ccb.dev.ea\_psi9000*), 104
- PSI9000Config (*class in hvl\_ccb.dev.ea\_psi9000*), 106
- PSI9000Error, 107
- PSI9000VisaCommunication (*class in hvl\_ccb.dev.ea\_psi9000*), 107
- PSI9000VisaCommunicationConfig (*class in hvl\_ccb.dev.ea\_psi9000*), 107
- PT100 (*LabJack.ThermocoupleType* attribute), 128
- PT1000 (*LabJack.ThermocoupleType* attribute), 128
- PT500 (*LabJack.ThermocoupleType* attribute), 128
- PTP (*ILS2T.Mode* attribute), 169
- PublicPropertiesReprMixin (*class in hvl\_ccb.dev.tiepie.utils*), 90
- PULSE (*TiePieGeneratorSignalType* attribute), 86
- pump\_init (*LaudaProRp245eConfig* attribute), 136
- PUMP\_LEVEL (*LaudaProRp245eCommand* attribute), 134
- ## Q
- QIL (*NewportConfigCommands* attribute), 142
- QUERY (*FuGProbusIVCommands* attribute), 114
- query() (*CryLasLaserSerialCommunication* method), 102
- query() (*FuGSerialCommunication* method), 118
- query() (*NewportSMC100PPSerialCommunication* method), 152
- query() (*PfeifferTPGSerialCommunication* method), 159
- query() (*SyncCommunicationProtocol* method), 14
- query() (*T560Communication* method), 54
- query() (*VisaCommunication* method), 30
- query\_all() (*CryLasLaserSerialCommunication* method), 102
- query\_command() (*LaudaProRp245eTcpCommunication* method), 137
- query\_multiple() (*NewportSMC100PPSerialCommunication* method), 152
- query\_polling() (*Luminox* method), 176
- query\_status() (*Technix* method), 78
- QUEUE (*AutoInstallMode* attribute), 53
- QUICK\_STOP (*SafetyStatus* attribute), 49
- QUICKSTOP (*ILS2T.State* attribute), 170
- QuickStop (*SafetyStatus* attribute), 70
- quickstop() (*ILS2T* method), 171
- quit (*Errors* attribute), 67
- quit\_error() (*BaseCube* method), 35
- quit\_error() (*Supercube2015Base* method), 60
- ## R
- R (*LabJack.ThermocoupleType* attribute), 128
- raise\_() (*FuGErrorcodes* method), 112
- RAMP\_ACC (*ILS2TRegAddr* attribute), 174
- RAMP\_CONTINUE (*LaudaProRp245eCommand* attribute), 134

- RAMP\_DECEL (*ILS2TRegAddr attribute*), 174
- RAMP\_DELETE (*LaudaProRp245eCommand attribute*), 134
- RAMP\_ITERATIONS (*LaudaProRp245eCommand attribute*), 134
- RAMP\_N\_MAX (*ILS2TRegAddr attribute*), 174
- RAMP\_PAUSE (*LaudaProRp245eCommand attribute*), 134
- RAMP\_SELECT (*LaudaProRp245eCommand attribute*), 134
- RAMP\_SET (*LaudaProRp245eCommand attribute*), 134
- RAMP\_START (*LaudaProRp245eCommand attribute*), 135
- RAMP\_STOP (*LaudaProRp245eCommand attribute*), 135
- RAMP\_TYPE (*ILS2TRegAddr attribute*), 174
- rampmode (*FuGProbusVSetRegisters property*), 117
- ramprate (*FuGProbusVSetRegisters property*), 117
- rampstate (*FuGProbusVSetRegisters property*), 117
- RAMPUPWARDS (*FuGRampModes attribute*), 117
- RATEDCURRENT (*FuGReadbackChannels attribute*), 117
- RATEDVOLTAGE (*FuGReadbackChannels attribute*), 117
- read() (*AsyncCommunicationProtocol method*), 11
- read() (*BaseCube method*), 35
- read() (*CryLasLaserSerialCommunication method*), 102
- read() (*LaudaProRp245eTcpCommunication method*), 137
- read() (*MBW973 method*), 139
- read() (*OpCuaCommunication method*), 21
- read() (*Supercube2015Base method*), 60
- read() (*Tcp method*), 26
- read\_all() (*AsyncCommunicationProtocol method*), 11
- read\_bytes() (*AsyncCommunicationProtocol method*), 12
- read\_bytes() (*SerialCommunication method*), 23
- read\_bytes() (*TelnetCommunication method*), 28
- read\_float() (*MBW973 method*), 139
- read\_holding\_registers() (*ModbusTcpCommunication method*), 18
- read\_input\_registers() (*ModbusTcpCommunication method*), 18
- read\_int() (*MBW973 method*), 139
- read\_measurement() (*RTO1024 method*), 164
- read\_measurements() (*MBW973 method*), 139
- read\_name() (*LJMCommunication method*), 15
- read\_nonempty() (*AsyncCommunicationProtocol method*), 12
- read\_output\_while\_polling (*TechnixConfig attribute*), 79
- read\_resistance() (*LabJack method*), 129
- read\_single\_bytes() (*SerialCommunication method*), 23
- read\_streaming() (*Luminos method*), 176
- read\_termination (*VisaCommunicationConfig attribute*), 32
- read\_text() (*AsyncCommunicationProtocol method*), 12
- read\_text() (*NewportSMC100PPSerialCommunication method*), 152
- read\_text\_nonempty() (*AsyncCommunicationProtocol method*), 12
- READ\_TEXT\_SKIP\_PREFIXES (*CryLasLaserSerialCommunication attribute*), 101
- read\_thermocouple() (*LabJack method*), 130
- readback\_data (*FuGProbusVConfigRegisters property*), 114
- READBACKCHANNEL (*FuGProbusIVCommands attribute*), 114
- ready (*BaseCube property*), 35
- READY (*CryLasLaser.AnswersStatus attribute*), 97
- READY (*ILS2T.State attribute*), 170
- READY (*NewportStates attribute*), 155
- ready() (*Supercube2015Base method*), 60
- READY\_ACTIVE (*CryLasLaser.LaserStatus attribute*), 98
- READY\_FROM\_DISABLE (*NewportSMC100PP.StateMessages attribute*), 144
- READY\_FROM\_HOMING (*NewportSMC100PP.StateMessages attribute*), 144
- READY\_FROM\_JOGGING (*NewportSMC100PP.StateMessages attribute*), 144
- READY\_FROM\_MOVING (*NewportSMC100PP.StateMessages attribute*), 144
- READY\_INACTIVE (*CryLasLaser.LaserStatus attribute*), 98
- record\_length (*TiePieOscilloscopeConfig property*), 90
- RED\_OPERATE (*SafetyStatus attribute*), 49
- RED\_READY (*SafetyStatus attribute*), 49
- RedOperate (*SafetyStatus attribute*), 71
- RedReady (*SafetyStatus attribute*), 71
- reg\_3 (*FuGProbusVDIRegisters property*), 115
- RegAddr (*ILS2T attribute*), 170
- RegDatatype (*ILS2T attribute*), 170
- register\_pulse\_time (*TechnixConfig attribute*), 79
- RELATIVE (*TiePieOscilloscopeTriggerLevelMode attribute*), 84
- RELATIVE\_POSITION\_MOTOR (*ILS2T.ActionsPtp attribute*), 169
- RELATIVE\_POSITION\_TARGET (*ILS2T.ActionsPtp attribute*), 169
- remote (*Technix property*), 78
- remove\_device() (*DeviceSequenceMixin method*), 93
- require\_block\_measurement\_support (*TiePieDeviceConfig attribute*), 81
- required\_keys() (*AsyncCommunicationProtocolCon-*

- fig* class method), 13
- required\_keys() (*BaseCubeConfiguration* class method), 37
- required\_keys() (*BaseCubeOpcUaCommunicationConfig* class method), 38
- required\_keys() (*CryLasAttenuatorConfig* class method), 96
- required\_keys() (*CryLasAttenuatorSerialCommunicationConfig* class method), 97
- required\_keys() (*CryLasLaserConfig* class method), 101
- required\_keys() (*CryLasLaserSerialCommunicationConfig* class method), 103
- required\_keys() (*EmptyConfig* class method), 94, 188
- required\_keys() (*FuGConfig* class method), 110
- required\_keys() (*FuGSerialCommunicationConfig* class method), 119
- required\_keys() (*HeinzingerConfig* class method), 120
- required\_keys() (*HeinzingerSerialCommunicationConfig* class method), 124
- required\_keys() (*ILS2TConfig* class method), 173
- required\_keys() (*ILS2TModbusTcpCommunicationConfig* class method), 174
- required\_keys() (*LaudaProRp245eConfig* class method), 136
- required\_keys() (*LaudaProRp245eTcpCommunicationConfig* class method), 138
- required\_keys() (*LJMCommunicationConfig* class method), 17
- required\_keys() (*LuminoxConfig* class method), 177
- required\_keys() (*LuminoxSerialCommunicationConfig* class method), 179
- required\_keys() (*MBW973Config* class method), 140
- required\_keys() (*MBW973SerialCommunicationConfig* class method), 141
- required\_keys() (*ModbusTcpCommunicationConfig* class method), 19
- required\_keys() (*NewportSMC100PPConfig* class method), 151
- required\_keys() (*NewportSMC100PPSerialCommunicationConfig* class method), 154
- required\_keys() (*OpcUaCommunicationConfig* class method), 22
- required\_keys() (*PfeifferTPGConfig* class method), 159
- required\_keys() (*PfeifferTPGSerialCommunicationConfig* class method), 160
- required\_keys() (*PICubeConfiguration* class method), 51
- required\_keys() (*PICubeOpcUaCommunicationConfig* class method), 52
- required\_keys() (*PSI9000Config* class method), 107
- required\_keys() (*PSI9000VisaCommunicationConfig* class method), 108
- required\_keys() (*RTO1024Config* class method), 168
- required\_keys() (*RTO1024VisaCommunicationConfig* class method), 168
- required\_keys() (*SerialCommunicationConfig* class method), 25
- required\_keys() (*SupercubeAOpcUaConfiguration* class method), 73
- required\_keys() (*SupercubeConfiguration* class method), 62
- required\_keys() (*SupercubeOpcUaCommunicationConfig* class method), 63
- required\_keys() (*T560CommunicationConfig* class method), 54
- required\_keys() (*T560Config* class method), 58
- required\_keys() (*TcpCommunicationConfig* class method), 27
- required\_keys() (*TechnixConfig* class method), 79
- required\_keys() (*TechnixSerialCommunicationConfig* class method), 75
- required\_keys() (*TechnixTelnetCommunicationConfig* class method), 76
- required\_keys() (*TelnetCommunicationConfig* class method), 29
- required\_keys() (*TiePieDeviceConfig* class method), 81
- required\_keys() (*VisaCommunicationConfig* class method), 32
- required\_keys() (*VisaDeviceConfig* class method), 182
- reset (*BreakdownDetection* attribute), 65
- RESET (*FuGProbusIVCommands* attribute), 114
- reset() (*FuGProbusIV* method), 113
- reset() (*NewportSMC100PP* method), 147
- reset() (*VisaDevice* method), 181
- reset\_error() (*ILS2T* method), 171
- reset\_interface() (*HeinzingerDI* method), 122
- reset\_ramp() (*LaudaProRp245e* method), 132
- resolution (*TiePieOscilloscopeConfig* property), 90
- response\_sleep\_time (*CryLasAttenuatorConfig* attribute), 96
- RISING (*TiePieOscilloscopeTriggerKind* attribute), 84
- RISING\_OR\_FALLING (*TiePieOscilloscopeTriggerKind* attribute), 84
- RMS\_CURRENT\_LIMIT (*NewportSMC100PP.MotorErrors* attribute), 143
- rpm\_max\_init (*ILS2TConfig* attribute), 173
- rs485\_address (*NewportSMC100PPConfig* attribute), 151
- RTO1024 (class in *hvl\_ccb.dev.rs\_rto1024*), 161
- RTO1024.TriggerModes (class in *hvl\_ccb.dev.rs\_rto1024*), 161

- RTO1024Config (class in *hvl\_ccb.dev.rs\_rto1024*), 167  
 RTO1024Error, 168  
 RTO1024VisaCommunication (class in *hvl\_ccb.dev.rs\_rto1024*), 168  
 RTO1024VisaCommunicationConfig (class in *hvl\_ccb.dev.rs\_rto1024*), 168  
 run() (*ExperimentManager* method), 190  
 run() (*LaudaProRp245e* method), 132  
 run\_continuous\_acquisition() (*RTO1024* method), 164  
 run\_single\_acquisition() (*RTO1024* method), 164  
 RUNNING (*ExperimentStatus* attribute), 190
- ## S
- S (*LabJack.ThermocoupleType* attribute), 128  
 SA (*NewportConfigCommands* attribute), 142  
 SafeGround (class in *hvl\_ccb.dev.tiepie.channel*), 82  
 Safety (class in *hvl\_ccb.dev.supercube2015.constants*), 70  
 SafetyStatus (class in *hvl\_ccb.dev.cube.constants*), 49  
 SafetyStatus (class in *hvl\_ccb.dev.supercube2015.constants*), 70  
 sample\_frequency (*TiePieOscilloscopeConfig* property), 90  
 save\_configuration() (*RTO1024* method), 164  
 save\_device\_configuration() (*T560* method), 57  
 save\_waveform\_history() (*RTO1024* method), 164  
 SCALE (*ILS2TRegAddr* attribute), 174  
 ScalingFactorValueError, 175  
 screw\_scaling (*NewportSMC100PPConfig* attribute), 151  
 send\_command() (*NewportSMC100PPSerialCommunication* method), 153  
 send\_command() (*PfeifferTPGSerialCommunication* method), 159  
 send\_hello() (*Client* method), 20  
 send\_pulses() (*LabJack* method), 130  
 send\_stop() (*NewportSMC100PPSerialCommunication* method), 153  
 Sensor (class in *hvl\_ccb.utils.conversion\_sensor*), 183  
 Sensor\_error (*PfeifferTPG.SensorStatus* attribute), 156  
 Sensor\_off (*PfeifferTPG.SensorStatus* attribute), 156  
 sensor\_status (*LuminoxMeasurementType* attribute), 177  
 SERIAL (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 135  
 serial\_number (*LuminoxMeasurementType* attribute), 177  
 serial\_number (*TiePieDeviceConfig* attribute), 81  
 SerialCommunication (class in *hvl\_ccb.comm.serial*), 23  
 SerialCommunicationBytesize (class in *hvl\_ccb.comm.serial*), 24  
 SerialCommunicationConfig (class in *hvl\_ccb.comm.serial*), 24  
 SerialCommunicationIOError, 25  
 SerialCommunicationParity (class in *hvl\_ccb.comm.serial*), 25  
 SerialCommunicationStopbits (class in *hvl\_ccb.comm.serial*), 25  
 Server (class in *hvl\_ccb.comm.opc*), 22  
 set\_acceleration() (*NewportSMC100PP* method), 147  
 set\_acquire\_length() (*RTO1024* method), 164  
 set\_ain\_differential() (*LabJack* method), 130  
 set\_ain\_range() (*LabJack* method), 130  
 set\_ain\_resistance() (*LabJack* method), 130  
 set\_ain\_resolution() (*LabJack* method), 130  
 set\_ain\_thermocouple() (*LabJack* method), 130  
 set\_analog\_output() (*LabJack* method), 131  
 set\_attenuation() (*CryLasAttenuator* method), 95  
 set\_cee16\_socket() (*Supercube2015Base* method), 60  
 set\_channel\_offset() (*RTO1024* method), 165  
 set\_channel\_position() (*RTO1024* method), 165  
 set\_channel\_range() (*RTO1024* method), 165  
 set\_channel\_scale() (*RTO1024* method), 165  
 set\_channel\_state() (*RTO1024* method), 166  
 set\_clock() (*LabJack* method), 131  
 set\_control\_mode() (*LaudaProRp245e* method), 132  
 set\_current() (*HeinzingerDI* method), 122  
 set\_current() (*HeinzingerPNC* method), 123  
 set\_digital\_output() (*LabJack* method), 131  
 set\_display\_unit() (*PfeifferTPG* method), 157  
 set\_earthing\_manual() (*Supercube2015Base* method), 60  
 set\_external\_temp() (*LaudaProRp245e* method), 133  
 set\_full\_scale\_mbar() (*PfeifferTPG* method), 157  
 set\_full\_scale\_unitless() (*PfeifferTPG* method), 158  
 set\_init\_attenuation() (*CryLasAttenuator* method), 95  
 set\_init\_shutter\_status() (*CryLasLaser* method), 99  
 set\_jog\_speed() (*ILS2T* method), 171  
 set\_lower\_limits() (*PSI9000* method), 105  
 set\_max\_acceleration() (*ILS2T* method), 171  
 set\_max\_deceleration() (*ILS2T* method), 171  
 set\_max\_rpm() (*ILS2T* method), 172  
 set\_measuring\_options() (*MBW973* method), 139  
 set\_message\_board() (*BaseCube* method), 35  
 set\_motor\_configuration() (*NewportSMC100PP* method), 147  
 set\_negative\_software\_limit() (*NewportSMC100PP* method), 148  
 set\_number\_of\_recordings() (*HeinzingerDI* method), 122

- set\_output() (*PSI9000 method*), 105  
 set\_positive\_software\_limit() (*NewportSMC100PP method*), 148  
 set\_pulse\_energy() (*CryLasLaser method*), 99  
 set\_pump\_level() (*LaudaProRp245e method*), 133  
 set\_ramp\_iterations() (*LaudaProRp245e method*), 133  
 set\_ramp\_program() (*LaudaProRp245e method*), 133  
 set\_ramp\_segment() (*LaudaProRp245e method*), 133  
 set\_ramp\_type() (*ILS2T method*), 172  
 set\_reference\_point() (*RTO1024 method*), 166  
 set\_register() (*FuGProbusV method*), 114  
 set\_remote\_control() (*Supercube2015Base method*), 60  
 set\_repetition\_rate() (*CryLasLaser method*), 99  
 set\_repetitions() (*RTO1024 method*), 166  
 set\_slope() (*Supercube2015WithFU method*), 72  
 set\_status\_board() (*BaseCube method*), 35  
 set\_support\_output() (*Supercube2015Base method*), 61  
 set\_support\_output\_impulse() (*Supercube2015Base method*), 61  
 set\_system\_lock() (*PSI9000 method*), 105  
 set\_t13\_socket() (*Supercube2015Base method*), 61  
 set\_target\_voltage() (*Supercube2015WithFU method*), 72  
 set\_temp\_set\_point() (*LaudaProRp245e method*), 133  
 set\_transmission() (*CryLasAttenuator method*), 95  
 set\_trigger\_level() (*RTO1024 method*), 166  
 set\_trigger\_mode() (*RTO1024 method*), 167  
 set\_trigger\_source() (*RTO1024 method*), 167  
 set\_upper\_limits() (*PSI9000 method*), 105  
 set\_voltage() (*HeinzingerDI method*), 122  
 set\_voltage() (*HeinzingerPNC method*), 123  
 set\_voltage\_current() (*PSI9000 method*), 105  
 SETCURRENT (*FuGProbusVRegisterGroups attribute*), 116  
 setup (*Power attribute*), 69  
 setvalue (*FuGProbusVSetRegisters property*), 117  
 SETVOLTAGE (*FuGProbusVRegisterGroups attribute*), 116  
 SEVENBITS (*SerialCommunicationBytesize attribute*), 24  
 SHORT\_CIRCUIT (*NewportSMC100PP.MotorErrors attribute*), 143  
 shunt (*LEM4000S attribute*), 183  
 SHUTDOWN\_CURRENT\_LIMIT (*PSI9000 attribute*), 104  
 SHUTDOWN\_VOLTAGE\_LIMIT (*PSI9000 attribute*), 104  
 ShutterStatus (*CryLasLaser attribute*), 98  
 ShutterStatus (*CryLasLaserConfig attribute*), 100  
 signal\_type (*TiePieGeneratorConfig property*), 86  
 SINE (*TiePieGeneratorSignalType attribute*), 86  
 SingleCommDevice (*class in hvl\_ccb.dev.base*), 94  
 SIXBITS (*SerialCommunicationBytesize attribute*), 24  
 SIXTEEN (*HeinzingerConfig.RecordingsEnum attribute*), 120  
 SIXTEEN\_BIT (*TiePieOscilloscopeResolution attribute*), 90  
 SL (*NewportConfigCommands attribute*), 142  
 SN (*FuGReadbackChannels attribute*), 117  
 SOFTWARE\_INTERNAL\_SIXTY (*CryLasLaser.RepetitionRates attribute*), 98  
 SOFTWARE\_INTERNAL\_TEN (*CryLasLaser.RepetitionRates attribute*), 98  
 SOFTWARE\_INTERNAL\_TWENTY (*CryLasLaser.RepetitionRates attribute*), 98  
 software\_revision (*LuminoxMeasurementType attribute*), 177  
 SPACE (*SerialCommunicationParity attribute*), 25  
 SPECIALRAMPUPWARDS (*FuGRampModes attribute*), 117  
 spoll() (*VisaCommunication method*), 31  
 spoll\_handler() (*VisaDevice method*), 181  
 SQUARE (*TiePieGeneratorSignalType attribute*), 86  
 SR (*NewportConfigCommands attribute*), 142  
 srq\_mask (*FuGProbusVConfigRegisters property*), 114  
 srq\_status (*FuGProbusVConfigRegisters property*), 115  
 stage\_configuration (*NewportSMC100PPConfig attribute*), 151  
 START (*LaudaProRp245eCommand attribute*), 135  
 start() (*BaseCube method*), 35  
 start() (*CryLasAttenuator method*), 95  
 start() (*CryLasLaser method*), 99  
 start() (*Device method*), 92  
 start() (*DeviceSequenceMixin method*), 93  
 start() (*ExperimentManager method*), 190  
 start() (*FuG method*), 109  
 start() (*FuGProbusIV method*), 113  
 start() (*HeinzingerDI method*), 122  
 start() (*HeinzingerPNC method*), 123  
 start() (*ILS2T method*), 172  
 start() (*LabJack method*), 131  
 start() (*LaudaProRp245e method*), 133  
 start() (*Luminox method*), 176  
 start() (*MBW973 method*), 139  
 start() (*NewportSMC100PP method*), 148  
 start() (*PfeifferTPG method*), 158  
 start() (*PSI9000 method*), 106  
 start() (*RTO1024 method*), 167  
 start() (*SingleCommDevice method*), 94  
 start() (*Supercube2015Base method*), 61  
 start() (*Technix method*), 78  
 start() (*TiePieGeneratorMixin method*), 86  
 start() (*TiePieI2CHostMixin method*), 87  
 start() (*TiePieOscilloscope method*), 89  
 start() (*VisaDevice method*), 181  
 start\_control() (*MBW973 method*), 139

- start\_measurement() (*TiePieOscilloscope method*), 89
- start\_polling() (*Poller method*), 180
- start\_ramp() (*LaudaProRp245e method*), 133
- STARTING (*ExperimentStatus attribute*), 190
- States (*NewportSMC100PP attribute*), 144
- status (*BaseCube property*), 35
- status (*ExperimentManager property*), 190
- status (*FuGProbusVConfigRegisters property*), 115
- status (*FuGProbusVDORegisters property*), 116
- status (*Technix property*), 78
- status\_1\_closed (*EarthingStick attribute*), 66
- status\_1\_connected (*EarthingStick attribute*), 66
- status\_1\_open (*EarthingStick attribute*), 66
- status\_2\_closed (*EarthingStick attribute*), 66
- status\_2\_connected (*EarthingStick attribute*), 66
- status\_2\_open (*EarthingStick attribute*), 66
- status\_3\_closed (*EarthingStick attribute*), 66
- status\_3\_connected (*EarthingStick attribute*), 66
- status\_3\_open (*EarthingStick attribute*), 66
- status\_4\_closed (*EarthingStick attribute*), 66
- status\_4\_connected (*EarthingStick attribute*), 66
- status\_4\_open (*EarthingStick attribute*), 66
- status\_5\_closed (*EarthingStick attribute*), 66
- status\_5\_connected (*EarthingStick attribute*), 66
- status\_5\_open (*EarthingStick attribute*), 66
- status\_6\_closed (*EarthingStick attribute*), 66
- status\_6\_connected (*EarthingStick attribute*), 66
- status\_6\_open (*EarthingStick attribute*), 66
- status\_closed() (*EarthingStick class method*), 66
- status\_connected() (*EarthingStick class method*), 66
- status\_error (*Safety attribute*), 70
- status\_green (*Safety attribute*), 70
- status\_open() (*EarthingStick class method*), 66
- status\_ready\_for\_red (*Safety attribute*), 70
- status\_red (*Safety attribute*), 70
- STATUSBYTE (*FuGReadbackChannels attribute*), 117
- stop (*Errors attribute*), 67
- STOP (*LaudaProRp245eCommand attribute*), 135
- stop() (*BaseCube method*), 35
- stop() (*CryLasLaser method*), 99
- stop() (*Device method*), 92
- stop() (*DeviceSequenceMixin method*), 93
- stop() (*ExperimentManager method*), 190
- stop() (*FuGProbusIV method*), 113
- stop() (*HeinzingerDI method*), 122
- stop() (*ILS2T method*), 172
- stop() (*LabJack method*), 131
- stop() (*LaudaProRp245e method*), 134
- stop() (*Luminox method*), 176
- stop() (*MBW973 method*), 139
- stop() (*NewportSMC100PP method*), 148
- stop() (*PfeifferTPG method*), 158
- stop() (*PSI9000 method*), 106
- stop() (*RTO1024 method*), 167
- stop() (*SingleCommDevice method*), 94
- stop() (*Supercube2015Base method*), 61
- stop() (*Technix method*), 78
- stop() (*TiePieGeneratorMixin method*), 86
- stop() (*TiePieI2CHostMixin method*), 87
- stop() (*TiePieOscilloscope method*), 89
- stop() (*VisaDevice method*), 181
- stop\_acquisition() (*RTO1024 method*), 167
- stop\_motion() (*NewportSMC100PP method*), 148
- stop\_number (*Errors attribute*), 67
- stop\_polling() (*Poller method*), 180
- stop\_ramp() (*LaudaProRp245e method*), 134
- STOP\_SAFETY\_STATUSES (*in module hvl\_ccb.dev.cube.constants*), 49
- stopbits (*CryLasAttenuatorSerialCommunicationConfig attribute*), 97
- stopbits (*CryLasLaserSerialCommunicationConfig attribute*), 103
- stopbits (*FuGSerialCommunicationConfig attribute*), 119
- stopbits (*HeinzingerSerialCommunicationConfig attribute*), 125
- stopbits (*LuminoxSerialCommunicationConfig attribute*), 179
- stopbits (*MBW973SerialCommunicationConfig attribute*), 141
- stopbits (*NewportSMC100PPSerialCommunicationConfig attribute*), 154
- stopbits (*PfeifferTPGSerialCommunicationConfig attribute*), 160
- Stopbits (*SerialCommunicationConfig attribute*), 24
- stopbits (*SerialCommunicationConfig attribute*), 25
- streaming (*LuminoxOutputMode attribute*), 178
- StrEnumBase (*class in hvl\_ccb.utils.enum*), 185
- sub\_handler (*BaseCubeOpcUaCommunicationConfig attribute*), 38
- sub\_handler (*OpcUaCommunicationConfig attribute*), 22
- sub\_handler (*SupercubeOpcUaCommunicationConfig attribute*), 63
- suitable\_range() (*TiePieOscilloscopeRange static method*), 83
- Supercube2015Base (*class in hvl\_ccb.dev.supercube2015.base*), 58
- Supercube2015WithFU (*class in hvl\_ccb.dev.supercube2015.typ\_a*), 71
- SupercubeAOpcUaCommunication (*class in hvl\_ccb.dev.supercube2015.typ\_a*), 72
- SupercubeAOpcUaConfiguration (*class in hvl\_ccb.dev.supercube2015.typ\_a*), 72
- SupercubeConfiguration (*class in hvl\_ccb.dev.supercube2015.base*), 61
- SupercubeOpcEndpoint (*class in*

- hvl\_ccb.dev.supercube2015.constants*), 71
- SupercubeOpcUaCommunication (class in *hvl\_ccb.dev.supercube2015.base*), 62
- SupercubeOpcUaCommunicationConfig (class in *hvl\_ccb.dev.supercube2015.base*), 62
- SupercubeSubscriptionHandler (class in *hvl\_ccb.dev.supercube2015.base*), 63
- switchto\_green (Safety attribute), 70
- switchto\_operate (Safety attribute), 70
- switchto\_ready (Safety attribute), 70
- SyncCommunicationProtocol (class in *hvl\_ccb.comm.base*), 14
- SyncCommunicationProtocolConfig (class in *hvl\_ccb.comm.base*), 14
- ## T
- T (*LabJack.ThermocoupleType* attribute), 128
- t13\_1 (*GeneralSockets* attribute), 67
- t13\_2 (*GeneralSockets* attribute), 67
- t13\_3 (*GeneralSockets* attribute), 67
- t13\_socket\_1 (*BaseCube* attribute), 36
- t13\_socket\_2 (*BaseCube* attribute), 36
- t13\_socket\_3 (*BaseCube* attribute), 36
- T13\_SOCKET\_PORTS (in module *hvl\_ccb.dev.supercube2015.constants*), 71
- T1MS (*FuGMonitorModes* attribute), 112
- T200MS (*FuGMonitorModes* attribute), 112
- T20MS (*FuGMonitorModes* attribute), 112
- T256US (*FuGMonitorModes* attribute), 112
- T4 (*LabJack.DeviceType* attribute), 127
- T4 (*LJMCommunicationConfig.DeviceType* attribute), 16
- T40MS (*FuGMonitorModes* attribute), 112
- T4MS (*FuGMonitorModes* attribute), 112
- T560 (class in *hvl\_ccb.dev.highland\_t560.device*), 56
- T560Communication (class in *hvl\_ccb.dev.highland\_t560.base*), 54
- T560CommunicationConfig (class in *hvl\_ccb.dev.highland\_t560.base*), 54
- T560Config (class in *hvl\_ccb.dev.highland\_t560.device*), 57
- T560Error, 55
- T7 (*LabJack.DeviceType* attribute), 127
- T7 (*LJMCommunicationConfig.DeviceType* attribute), 16
- T7\_PRO (*LabJack.DeviceType* attribute), 127
- T7\_PRO (*LJMCommunicationConfig.DeviceType* attribute), 16
- T800MS (*FuGMonitorModes* attribute), 112
- T80MS (*FuGMonitorModes* attribute), 113
- target\_pulse\_energy (*CryLasLaser* property), 100
- Tcp (class in *hvl\_ccb.comm.tcp*), 26
- TCP (*LJMCommunicationConfig.ConnectionType* attribute), 16
- TcpCommunicationConfig (class in *hvl\_ccb.comm.tcp*), 27
- TCPIP\_INSTR (*VisaCommunicationConfig.InterfaceType* attribute), 31
- TCPIP\_SOCKET (*VisaCommunicationConfig.InterfaceType* attribute), 31
- TEC1 (*CryLasLaser.AnswersStatus* attribute), 97
- TEC2 (*CryLasLaser.AnswersStatus* attribute), 98
- Technix (class in *hvl\_ccb.dev.technix.device*), 77
- TechnixConfig (class in *hvl\_ccb.dev.technix.device*), 78
- TechnixError, 74
- TechnixFaultError, 74
- TechnixSerialCommunication (class in *hvl\_ccb.dev.technix.base*), 74
- TechnixSerialCommunicationConfig (class in *hvl\_ccb.dev.technix.base*), 74
- TechnixTelnetCommunication (class in *hvl\_ccb.dev.technix.base*), 75
- TechnixTelnetCommunicationConfig (class in *hvl\_ccb.dev.technix.base*), 75
- TelnetCommunication (class in *hvl\_ccb.comm.telnet*), 28
- TelnetCommunicationConfig (class in *hvl\_ccb.comm.telnet*), 28
- TelnetError, 29
- TEMP (*ILS2TRegAddr* attribute), 174
- TEMP\_SET\_POINT (*LaudaProRp245eCommand* attribute), 135
- temp\_set\_point\_init (*LaudaProRp245eConfig* attribute), 137
- Temperature (class in *hvl\_ccb.utils.conversion\_unit*), 184
- temperature\_sensor (*LuminoxMeasurementType* attribute), 178
- temperature\_unit (*LMT70A* attribute), 183
- TEN (*LabJack.AInRange* attribute), 126
- TEN (*LabJack.CalMicroAmpere* attribute), 126
- TEN\_MHZ (*LabJack.ClockFrequency* attribute), 127
- terminator (*AsyncCommunicationProtocolConfig* attribute), 14
- terminator (*CryLasAttenuatorSerialCommunicationConfig* attribute), 97
- terminator (*CryLasLaserSerialCommunicationConfig* attribute), 103
- TERMINATOR (*FuGProbusIVCommands* attribute), 114
- terminator (*FuGProbusVConfigRegisters* property), 115
- terminator (*FuGSerialCommunicationConfig* attribute), 119
- terminator (*HeinzingerSerialCommunicationConfig* attribute), 125
- terminator (*LaudaProRp245eTcpCommunicationConfig* attribute), 138
- terminator (*LuminoxSerialCommunicationConfig* attribute), 179
- terminator (*MBW973SerialCommunicationConfig* at-

tribute), 141

terminator (*NewportSMC100PPSerialCommunicationConfig* attribute), 154

terminator (*PfeifferTPGSerialCommunicationConfig* attribute), 160

terminator (*T560CommunicationConfig* attribute), 55

terminator\_str() (*SerialCommunicationConfig* method), 25

THIRTY\_TWO\_BIT (*LabJack.BitLimit* attribute), 126

TiePieDeviceConfig (class in *hvl\_ccb.dev.tiepie.base*), 80

TiePieDeviceType (class in *hvl\_ccb.dev.tiepie.base*), 81

TiePieError, 81

TiePieGeneratorConfig (class in *hvl\_ccb.dev.tiepie.generator*), 85

TiePieGeneratorConfigLimits (class in *hvl\_ccb.dev.tiepie.generator*), 86

TiePieGeneratorMixin (class in *hvl\_ccb.dev.tiepie.generator*), 86

TiePieGeneratorSignalType (class in *hvl\_ccb.dev.tiepie.generator*), 86

TiePieHS5 (class in *hvl\_ccb.dev.tiepie.device*), 84

TiePieHS6 (class in *hvl\_ccb.dev.tiepie.device*), 85

TiePieI2CHostConfig (class in *hvl\_ccb.dev.tiepie.i2c*), 87

TiePieI2CHostConfigLimits (class in *hvl\_ccb.dev.tiepie.i2c*), 87

TiePieI2CHostMixin (class in *hvl\_ccb.dev.tiepie.i2c*), 87

TiePieOscilloscope (class in *hvl\_ccb.dev.tiepie.oscilloscope*), 88

TiePieOscilloscopeAutoResolutionModes (class in *hvl\_ccb.dev.tiepie.oscilloscope*), 89

TiePieOscilloscopeChannelConfig (class in *hvl\_ccb.dev.tiepie.channel*), 82

TiePieOscilloscopeChannelConfigLimits (class in *hvl\_ccb.dev.tiepie.channel*), 83

TiePieOscilloscopeChannelCoupling (class in *hvl\_ccb.dev.tiepie.channel*), 83

TiePieOscilloscopeConfig (class in *hvl\_ccb.dev.tiepie.oscilloscope*), 89

TiePieOscilloscopeConfigLimits (class in *hvl\_ccb.dev.tiepie.oscilloscope*), 90

TiePieOscilloscopeRange (class in *hvl\_ccb.dev.tiepie.channel*), 83

TiePieOscilloscopeResolution (class in *hvl\_ccb.dev.tiepie.oscilloscope*), 90

TiePieOscilloscopeTriggerKind (class in *hvl\_ccb.dev.tiepie.channel*), 83

TiePieOscilloscopeTriggerLevelMode (class in *hvl\_ccb.dev.tiepie.channel*), 84

TiePieWS5 (class in *hvl\_ccb.dev.tiepie.device*), 85

timeout (*CryLasAttenuatorSerialCommunicationConfig* attribute), 97

timeout (*CryLasLaserSerialCommunicationConfig* attribute), 103

timeout (*FuGSerialCommunicationConfig* attribute), 119

timeout (*HeinzingerSerialCommunicationConfig* attribute), 125

timeout (*LuminosSerialCommunicationConfig* attribute), 179

timeout (*MBW973SerialCommunicationConfig* attribute), 141

timeout (*NewportSMC100PPSerialCommunicationConfig* attribute), 154

timeout (*PfeifferTPGSerialCommunicationConfig* attribute), 160

timeout (*SerialCommunicationConfig* attribute), 25

timeout (*TelnetCommunicationConfig* attribute), 29

timeout (*VisaCommunicationConfig* attribute), 32

timeout\_interval (*BaseCubeConfiguration* attribute), 37

timeout\_status\_change (*BaseCubeConfiguration* attribute), 37

timeout\_test\_parameters (*PICubeConfiguration* attribute), 51

Torr (*PfeifferTPG.PressureUnits* attribute), 155

TORR (*Pressure* attribute), 184

TPG25xA (*PfeifferTPGConfig.Model* attribute), 158

TPGx6x (*PfeifferTPGConfig.Model* attribute), 158

TPR (*PfeifferTPG.SensorTypes* attribute), 156

transmission (*CryLasAttenuator* property), 95

TRIANGLE (*TiePieGeneratorSignalType* attribute), 86

trigger\_enabled (*TiePieOscilloscopeChannelConfig* property), 83

trigger\_hysteresis (*TiePieOscilloscopeChannelConfig* property), 83

trigger\_kind (*TiePieOscilloscopeChannelConfig* property), 83

trigger\_level (*T560* property), 57

trigger\_level (*TiePieOscilloscopeChannelConfig* property), 83

trigger\_level\_mode (*TiePieOscilloscopeChannelConfig* property), 83

trigger\_mode (*T560* property), 57

trigger\_timeout (*TiePieOscilloscopeConfig* property), 90

triggered (*BreakdownDetection* attribute), 65

TriggerMode (class in *hvl\_ccb.dev.highland\_t560.base*), 55

TWELVE\_BIT (*TiePieOscilloscopeResolution* attribute), 90

TWELVE\_HUNDRED\_FIFTY\_KHZ (*LabJack.ClockFrequency* attribute), 127

TWENTY\_FIVE\_HUNDRED\_KHZ (*LabJack.ClockFrequency* attribute), 127

TWENTY\_MHZ (*LabJack.ClockFrequency* attribute), 127  
 TWENTY\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
 TWO (*HeinzingerConfig.RecordingsEnum* attribute), 120  
 TWO (*SerialCommunicationStopbits* attribute), 26  
 TWO\_HUNDRED (*LabJack.CalMicroAmpere* attribute), 126  
 TWO\_HUNDRED\_MILLI\_VOLT (*TiePieOscilloscopeRange* attribute), 83  
 TWO\_VOLT (*TiePieOscilloscopeRange* attribute), 83

## U

Underrange (*PfeifferTPG.SensorStatus* attribute), 156  
 Unit (*class in hvl\_ccb.utils.conversion\_unit*), 184  
 unit (*ILS2TModbusTcpCommunicationConfig* attribute), 174  
 unit (*ModbusTcpCommunicationConfig* attribute), 19  
 unit (*PfeifferTPG* property), 158  
 unit\_current (*HeinzingerPNC* property), 123  
 unit\_voltage (*HeinzingerPNC* property), 123  
 UNKNOWN (*HeinzingerDI.OutputStatus* attribute), 121  
 UNKNOWN (*HeinzingerPNC.UnitCurrent* attribute), 122  
 UNKNOWN (*HeinzingerPNC.UnitVoltage* attribute), 123  
 UNKNOWN (*TiePieGeneratorSignalType* attribute), 86  
 UNKNOWN (*TiePieOscilloscopeAutoResolutionModes* attribute), 89  
 UNKNOWN (*TiePieOscilloscopeTriggerLevelMode* attribute), 84  
 UNREADY\_INACTIVE (*CryLasLaser.LaserStatus* attribute), 98  
 update\_laser\_status() (*CryLasLaser* method), 100  
 update\_parameter (*OpCuaCommunicationConfig* attribute), 22  
 update\_repetition\_rate() (*CryLasLaser* method), 100  
 update\_shutter\_status() (*CryLasLaser* method), 100  
 update\_target\_pulse\_energy() (*CryLasLaser* method), 100  
 UpdateEspStageInfo (*NewportSMC100PPConfig.EspStageConfig* attribute), 149  
 UPPER\_TEMP (*LaudaProRp245eCommand* attribute), 135  
 upper\_temp (*LaudaProRp245eConfig* attribute), 137  
 USB (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 135  
 USB (*LJMCommunicationConfig.ConnectionType* attribute), 16  
 use\_external\_clock() (*T560* method), 57  
 user\_position\_offset (*NewportSMC100PPConfig* attribute), 151  
 user\_steps() (*ILS2T* method), 172

## V

V (*HeinzingerPNC.UnitVoltage* attribute), 123  
 VA (*NewportConfigCommands* attribute), 142

validate\_and\_resolve\_host() (*in module hvl\_ccb.utils.validation*), 186  
 validate\_bool() (*in module hvl\_ccb.utils.validation*), 186  
 validate\_number() (*in module hvl\_ccb.utils.validation*), 186  
 validate\_pump\_level() (*LaudaProRp245e* method), 134  
 validate\_tcp\_port() (*in module hvl\_ccb.utils.validation*), 187  
 value (*FuGProbusVMonitorRegisters* property), 116  
 value (*LabJack.AInRange* property), 126  
 value\_raw (*FuGProbusVMonitorRegisters* property), 116  
 ValueEnum (*class in hvl\_ccb.utils.enum*), 185  
 VB (*NewportConfigCommands* attribute), 142  
 velocity (*NewportSMC100PPConfig* attribute), 151  
 visa\_backend (*VisaCommunicationConfig* attribute), 32  
 VisaCommunication (*class in hvl\_ccb.comm.visa*), 30  
 VisaCommunicationConfig (*class in hvl\_ccb.comm.visa*), 31  
 VisaCommunicationConfig.InterfaceType (*class in hvl\_ccb.comm.visa*), 31  
 VisaCommunicationError, 32  
 VisaDevice (*class in hvl\_ccb.dev.visa*), 180  
 VisaDeviceConfig (*class in hvl\_ccb.dev.visa*), 181  
 VOLT (*ILS2TRegAddr* attribute), 174  
 Volt (*PfeifferTPG.PressureUnits* attribute), 155  
 voltage (*FuG* property), 110  
 VOLTAGE (*FuGProbusIVCommands* attribute), 114  
 VOLTAGE (*FuGReadbackChannels* attribute), 117  
 voltage (*Technix* property), 78  
 voltage\_actual (*PICube* property), 51  
 voltage\_lower\_limit (*PSI9000Config* attribute), 107  
 voltage\_max (*PICube* property), 51  
 voltage\_max (*Power* attribute), 69  
 voltage\_monitor (*FuG* property), 110  
 voltage\_primary (*PICube* property), 51  
 voltage\_primary (*Power* attribute), 69  
 voltage\_regulation (*Technix* property), 78  
 voltage\_slope (*Power* attribute), 69  
 voltage\_target (*Power* attribute), 69  
 voltage\_upper\_limit (*PSI9000Config* attribute), 107

## W

WAIT\_AFTER\_WRITE (*VisaCommunication* attribute), 30  
 wait\_for\_polling\_result() (*Poller* method), 180  
 wait\_operation\_complete() (*VisaDevice* method), 181  
 wait\_sec\_initialisation (*PSI9000Config* attribute), 107  
 wait\_sec\_max\_disable (*ILS2TConfig* attribute), 173  
 wait\_sec\_post\_absolute\_position (*ILS2TConfig* attribute), 173

`wait_sec_post_activate` (*LuminosConfig* attribute), 177

`wait_sec_post_cannot_disable` (*ILS2TConfig* attribute), 173

`wait_sec_post_enable` (*ILS2TConfig* attribute), 173

`wait_sec_post_relative_step` (*ILS2TConfig* attribute), 173

`wait_sec_pre_read_or_write` (*LaudaProRp245eTcpCommunicationConfig* attribute), 138

`wait_sec_read_text_nonempty` (*AsyncCommunicationProtocolConfig* attribute), 14

`wait_sec_read_text_nonempty` (*FuGSerialCommunicationConfig* attribute), 119

`wait_sec_read_text_nonempty` (*HeinzingerSerialCommunicationConfig* attribute), 125

`wait_sec_retry_get_device` (*TiePieDeviceConfig* attribute), 81

`wait_sec_settings_effect` (*PSI9000Config* attribute), 107

`wait_sec_stop_commands` (*FuGConfig* attribute), 110

`wait_sec_stop_commands` (*HeinzingerConfig* attribute), 120

`wait_sec_system_lock` (*PSI9000Config* attribute), 107

`wait_sec_trials_activate` (*LuminosConfig* attribute), 177

`wait_timeout_retry_sec` (*OpcUaCommunicationConfig* attribute), 22

`wait_until_motor_initialized()` (*NewportSMC100PP* method), 149

`wait_until_ready()` (*CryLasLaser* method), 100

`waveform` (*TiePieGeneratorConfig* property), 86

`WIFI` (*LJMCommunicationConfig.ConnectionType* attribute), 16

`wrap_libtiepie_exception()` (in module *hvl\_ccb.dev.tiepie.base*), 81

`write()` (*AsyncCommunicationProtocol* method), 12

`write()` (*BaseCube* method), 36

`write()` (*MBW973* method), 139

`write()` (*OpcUaCommunication* method), 21

`write()` (*Supercube2015Base* method), 61

`write()` (*Tcp* method), 26

`write()` (*VisaCommunication* method), 31

`write_absolute_position()` (*ILS2T* method), 172

`write_bytes()` (*AsyncCommunicationProtocol* method), 12

`write_bytes()` (*SerialCommunication* method), 23

`write_bytes()` (*TelnetCommunication* method), 28

`write_command()` (*LaudaProRp245eTcpCommunication* method), 137

`write_name()` (*LJMCommunication* method), 16

`write_names()` (*LJMCommunication* method), 16

`write_registers()` (*ModbusTcpCommunication* method), 18

`write_relative_step()` (*ILS2T* method), 172

`write_termination` (*VisaCommunicationConfig* attribute), 32

`write_text()` (*AsyncCommunicationProtocol* method), 13

`WRONG_ESP_STAGE` (*NewportSMC100PP.MotorErrors* attribute), 143

## X

`x_stat` (*FuGProbusVDIRegisters* property), 115

`XOUTPUTS` (*FuGProbusIVCommands* attribute), 114

## Y

`YES` (*FuGDigitalVal* attribute), 111

## Z

`ZX` (*NewportConfigCommands* attribute), 142