
HVL Common Code Base Documentation

Release 0.14.4

Henning Janssen, Chi-Ching Hsu

Dec 22, 2023

CONTENTS:

1	HVL Common Code Base	1
1.1	Features	1
1.2	Documentation	5
1.3	Credits	5
2	Installation	7
2.1	Stable release	7
2.2	From sources	7
2.3	Additional system libraries	8
3	Usage	9
4	API Documentation	11
4.1	hvl_ccb	11
5	Contributing	205
5.1	Types of Contributions	205
5.2	Get Started!	206
5.3	Merge Request Guidelines	207
5.4	Tips	207
5.5	Deploying	208
6	Credits	211
6.1	Active Maintainers	211
6.2	Authors and Contributors	211
7	History	213
7.1	0.14.4 (2023-12-22)	213
7.2	0.14.3 (2023-11-17)	213
7.3	0.14.2 (2023-09-07)	213
7.4	0.14.1 (2023-08-21)	214
7.5	0.14.0 (2023-07-28)	214
7.6	0.13.3 (2023-03-31)	214
7.7	0.13.2 (2023-03-17)	215
7.8	0.13.1 (2023-03-03)	215
7.9	0.13.0 (2023-01-27)	215
7.10	0.12.3 (2022-12-27)	215
7.11	0.12.2 (2022-11-29)	216
7.12	0.12.1 (2022-10-31)	216
7.13	0.12.0 (2022-10-17)	216
7.14	0.11.1 (2022-09-15)	216

7.15	0.11.0 (2022-06-22)	217
7.16	0.10.3 (2022-03-21)	217
7.17	0.10.2 (2022-02-28)	217
7.18	0.10.1 (2022-01-24)	217
7.19	0.10.0 (2022-01-17)	218
7.20	0.9.0 (2022-01-07)	218
7.21	0.8.5 (2021-11-05)	219
7.22	0.8.4 (2021-10-22)	219
7.23	0.8.3 (2021-09-27)	219
7.24	0.8.2 (2021-08-27)	219
7.25	0.8.1 (2021-08-13)	220
7.26	0.8.0 (2021-07-02)	220
7.27	0.7.1 (2021-06-04)	220
7.28	0.7.0 (2021-05-25)	220
7.29	0.6.1 (2021-05-08)	221
7.30	0.6.0 (2021-04-23)	221
7.31	0.5.0 (2020-11-11)	222
7.32	0.4.0 (2020-07-16)	222
7.33	0.3.5 (2020-02-18)	223
7.34	0.3.4 (2019-12-20)	223
7.35	0.3.3 (2019-05-08)	223
7.36	0.3.2 (2019-05-08)	223
7.37	0.3.1 (2019-05-02)	224
7.38	0.3 (2019-05-02)	224
7.39	0.2.1 (2019-04-01)	224
7.40	0.2.0 (2019-03-31)	224
7.41	0.1.0 (2019-02-06)	224
8	Indices and tables	225
	Python Module Index	227
	Index	229

**CHAPTER
ONE**

HVL COMMON CODE BASE

Python common code base (CCB) to control devices, which are used in high-voltage research. All implemented devices are used and tested in the High Voltage Laboratory ([HVL](#)) of the Federal Institute of Technology Zurich (ETH Zurich).

- Free software: GNU General Public License v3
- Copyright (c) 2019-2023 ETH Zurich, SIS ID and HVL D-ITET

1.1 Features

For managing multi-device experiments instantiate the `ExperimentManager` utility class.

1.1.1 Devices

The device wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, various settings and options, as well as start/stop methods. Currently wrappers are available to control the following devices:

Function/Type	Devices
Bench Multimeter	Fluke 8845A and 8846A 6.5 Digit Precision Multimeter
Data acquisition	LabJack (T4, T7, T7-PRO; requires LJM Library) Pico Technology PT-104 Platinum Resistance Data Logger (requires PicoSDK/libusbpt104)
Digital Delay Generator	Highland T560
Digital IO	LabJack (T4, T7, T7-PRO; requires LJM Library)
Experiment control	HVL Cube with and without Power Inverter
Gas Analyser	MBW 973-SF6 gas dew point mirror analyzer Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges SST Luminox oxygen sensor
Laser	CryLaS pulsed laser CryLaS laser attenuator
Oscilloscope	Rhode & Schwarz RTO 1024 TiePie (HS5, HS6, WS5)
Power supply	Elektro-Automatik PSI9000 FuG Elektronik Heinzinger PNC Technix capacitor charger
Stepper motor drive	Newport SMC100PP Schneider Electric ILS2T
Temperature control	Lauda PRO RP 245 E circulation thermostat
Waveform generator	Chapter 1. HVL Common Code Base TiePie (HS5, WS5)

Each device uses at least one standardised communication protocol wrapper.

1.1.2 Communication protocols

In hvl_ccb by “communication protocol” we mean different levels of communication standards, from the low level actual communication protocols like serial communication to application level interfaces like VISA TCP standard. There are also devices in hvl_ccb that use a dummy communication protocol; this is because these devices are build on proprietary manufacturer libraries that communicate with the corresponding devices, as in the case of TiePie or LabJack devices.

The communication protocol wrappers in hvl_ccb provide a standardised API with configuration dataclasses, as well as open/close and read/write/query methods. Currently, wrappers for the following communication protocols are available:

Communication protocol	Devices using
Modbus TCP	Schneider Electric ILS2T stepper motor drive
OPC UA	HVL Cube with and without Power Inverter
Serial	CryLaS pulsed laser and laser attenuator FuG Elektronik power supply (e.g. capacitor charger HCK) using the Probus V protocol Heinzinger PNC power supply using Heinzinger Digital Interface I/II SST Luminox oxygen sensor MBW 973-SF6 gas dew point mirror analyzer Newport SMC100PP single axis driver for 2-phase stepper motors Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges Technix capacitor charger
TCP	Lauda PRO RP 245 E circulation thermostat
Telnet	Technix capacitor charger Fluke 8845A and 8846
VISA TCP	Elektro-Automatik PSI9000 DC power supply Rhode & Schwarz RTO 1024 oscilloscope
<i>proprietary</i>	LabJack (T4, T7, T7-PRO) devices, which communicate via LJM Library Pico Technology PT-104 Platinum Resistance Data Logger, which communicate via PicoSDK/libusbpt104 TiePie (HS5, HS6, WS5) oscilloscopes and generators, which communicate via LibTiePie SDK

1.1.3 Sensor and Unit Conversion Utility

The Conversion Utility is a submodule that allows on the one hand a unified implementation of hardware-sensors and on the other hand provides a unified way to convert units. Furthermore it is possible to map two ranges on to each other. This can be useful to convert between for example and 4 - 20 mA and 0 - 10 V, both of them are common as sensor out- or input. Moreover, a subclass allows the mapping of a bit-range to any other range. For example a 12 bit number (0-4095) to 0 - 10. All utilities can be used with single numbers (`int`, `float`) as well as array-like structures containing single numbers (`np.array()`, `list`, `dict`, `tuple`).

Currently the following sensors are implemented:

- LEM LT 4000S
- LMT 70A

The following unit conversion classes are implemented:

- Temperature (Kelvin, Celsius, Fahrenheit)
- Pressure (Pascal, Bar, Atmosphere, Psi, Torr, Millimeter Mercury)

1.2 Documentation

Note: if you're planning to contribute to the `hvl_ccb` project read the **Contributing** section in the HVL CCB documentation.

Do either:

- read [HVL CCB documentation at RTD](#),

or

- build and read HVL CCB documentation locally; install first `Graphviz` (make sure to have the `dot` command in the executable search path) and the Python build requirements for documentation:

```
$ pip install docs/requirements.txt
```

and then either on Windows in Git BASH run:

```
$ ./make.sh docs
```

or from any other shell with GNU Make installed run:

```
$ make docs
```

The target index HTML ("`docs/_build/html/index.html`") should open automatically in your Web browser.

1.3 Credits

This package was created with [Cookiecutter](#) and the [audreyr/cookiecutter-pypackage](#) project template.

INSTALLATION

2.1 Stable release

To install HVL Common Code Base, run this command in your terminal:

```
$ pip install hvl_ccb
```

To install HVL Common Code Base with optional Python libraries that require manual installations of additional system libraries, you need to specify on installation extra requirements corresponding to these controllers. For instance, to install Python requirements for LabJack and TiePie devices, run:

```
$ pip install "hvl_ccb[tiepie,labjack]"
```

See below for the info about additional system libraries and the corresponding extra requirements.

To install all extra requirements run:

```
$ pip install "hvl_ccb[all]"
```

This is the preferred method to install HVL Common Code Base, as it will always install the most recent stable release.

If you don't have `pip` installed, this [Python installation guide](#) can guide you through the process.

2.2 From sources

The sources for HVL Common Code Base can be downloaded from the [GitLab](#) repo.

You can either clone the repository:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

Or download the [tarball](#):

```
$ curl -OL https://gitlab.com/ethz_hvl/hvl_ccb/-/archive/master/hvl_ccb.tar.gz
```

Once you have a copy of the source, you can install it with:

```
$ pip install .
```

2.3 Additional system libraries

If you have installed `hvl_ccb` with any of the extra features corresponding to device controllers, you must additionally install respective system library; these are:

Extra feature	Additional system library
<code>labjack</code>	LJM Library
<code>picotech</code>	PicoSDK (Windows) / libusbpt104 (Ubuntu/Debian)

For more details on installation of the libraries see docstrings of the corresponding `hvl_ccb` modules.

**CHAPTER
THREE**

USAGE

To use HVL Common Code Base in a project:

```
import hvl_ccb
```


API DOCUMENTATION


4.1 hvl_ccb

4.1.1 Subpackages

[hvl_ccb.comm](#)

Submodules

[hvl_ccb.comm.base](#)

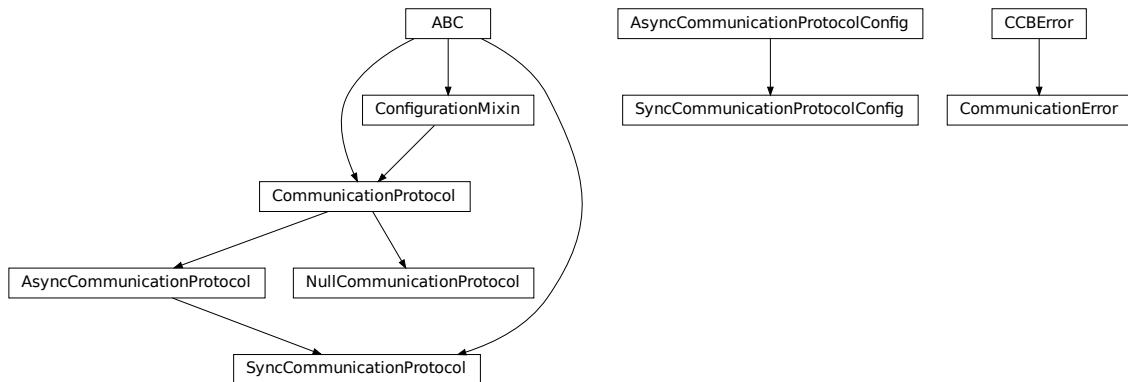

Module with base classes for communication protocols.

```
class AsyncCommunicationProtocol(config)
    Bases: CommunicationProtocol
    Abstract base class for asynchronous communication protocols
    static config_cls() → type[AsyncCommunicationProtocolConfig]
        Return the default configdataclass class.
```

Returns

a reference to the default configdataclass class

read() → str

Read a single line of text as *str* from the communication.

Returns

text as *str* including the terminator, which can also be empty “”

read_all(*n_attempts_max*: int | None = None, *attempt_interval_sec*: int | float | None = None) → str | *None*

Read all lines of text from the connection till nothing is left to read.

Parameters

- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt_interval_sec** – time between the reading attempts

Returns

A multi-line *str* including the terminator internally

abstract read_bytes() → bytes

Read a single line as *bytes* from the communication.

This method uses *self.access_lock* to ensure thread-safety.

Returns

a single line as *bytes* containing the terminator, which can also be empty b””

read_nonempty(*n_attempts_max*: int | None = None, *attempt_interval_sec*: int | float | None = None) → str | *None*

Try to read a non-empty single line of text as *str* from the communication. If the host does not reply or reply with white space only, it will return None.

Returns

a non-empty text as a *str* or *None* in case of an empty string

Parameters

- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt_interval_sec** – time between the reading attempts

read_text() → str

Read one line of text from the serial port. The input buffer may hold additional data afterwards, since only one line is read.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Returns

String read from the serial port; “ if there was nothing to read.

Raises

SerialCommunicationIOError – when communication port is not opened

read_text_nonempty(*n_attempts_max*: int | None = None, *attempt_interval_sec*: int | float | None = None)
→ str | *None*

Reads from the serial port, until a non-empty line is found, or the number of attempts is exceeded.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Attention: in contrast to *read_text*, the returned answer will be stripped of a whitespace newline terminator at the end, if such terminator is set in the initial configuration (default).

Parameters

- **n_attempts_max** – maximum number of read attempts

- **attempt_interval_sec** – time between the reading attempts

Returns

String read from the serial port; ‘’ if number of attempts is exceeded or serial port is not opened.

write(*text: str*)

Write text as *str* to the communication.

Parameters

text – test as a *str* to be written

abstract write_bytes(*data: bytes*) → int

Write data as *bytes* to the communication.

This method uses *self.access_lock* to ensure thread-safety.

Parameters

data – data as *bytes*-string to be written

Returns

number of bytes written

write_text(*text: str*)

Write text to the serial port. The text is encoded and terminated by the configured terminator.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Parameters

text – Text to send to the port.

Raises

SerialCommunicationIOError – when communication port is not opened

```
class AsyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: int | float = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10)
```

Bases: *object*

Base configuration data class for asynchronous communication protocols

clean_values()**default_n_attempts_read_text_nonempty: int = 10**

default number of attempts to read a non-empty text

encoding: str = 'utf-8'

Standard encoding of the connection. Typically this is *utf-8*, but can also be *latin-1* or something from here: <https://docs.python.org/3/library/codecs.html#standard-encodings>

encoding_error_handling: str = 'strict'

Encoding error handling scheme as defined here: <https://docs.python.org/3/library/codecs.html#error-handlers> By default strict error handling that raises *UnicodeError*.

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

terminator: bytes = b'\r\n'

The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations. This defines the end of a single line.

wait_sec_read_text_nonempty: int | float = 0.5

time to wait between attempts of reading a non-empty text

exception CommunicationError

Bases: *CCBError*

class CommunicationProtocol(config)

Bases: *ConfigurationMixin*, ABC

Communication protocol abstract base class.

Specifies the methods to implement for communication protocol, as well as implements some default settings and checks.

access_lock

Access lock to use with context manager when accessing the communication protocol (thread safety)

abstract close()

Close the communication protocol

abstract open()

Open communication protocol

class NullCommunicationProtocol(config)

Bases: *CommunicationProtocol*

Communication protocol that does nothing.

close() → None

Void close function.

static config_cls() → type[*EmptyConfig*]

Empty configuration

Returns

EmptyConfig

open() → *None*

Void open function.

class SyncCommunicationProtocol(*config*)

Bases: *AsyncCommunicationProtocol*, ABC

Abstract base class for synchronous communication protocols with *query()*

static config_cls() → type[*SyncCommunicationProtocolConfig*]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

query(*command*: str, *n_attempts_max*: int | *None* = *None*, *attempt_interval_sec*: int | float | *None* = *None*)

→ str | *None*

Send a command to the interface and handle the status message. Possibly raises an exception.

Parameters

- **command** – Command to send
- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read as answer
- **attempt_interval_sec** – time between the reading attempts

Returns

Answer from the interface, which can be None instead of an empty reply

class SyncCommunicationProtocolConfig(*terminator*: bytes = b'\r\n', *encoding*: str = 'utf-8',

encoding_error_handling: str = 'strict',


wait_sec_read_text_nonempty: int | float = 0.5,

default_n_attempts_read_text_nonempty: int = 10)

Bases: *AsyncCommunicationProtocolConfig*

Base configuration data class for synchronous communication protocols

hvl_ccb.comm.labjack_ljm

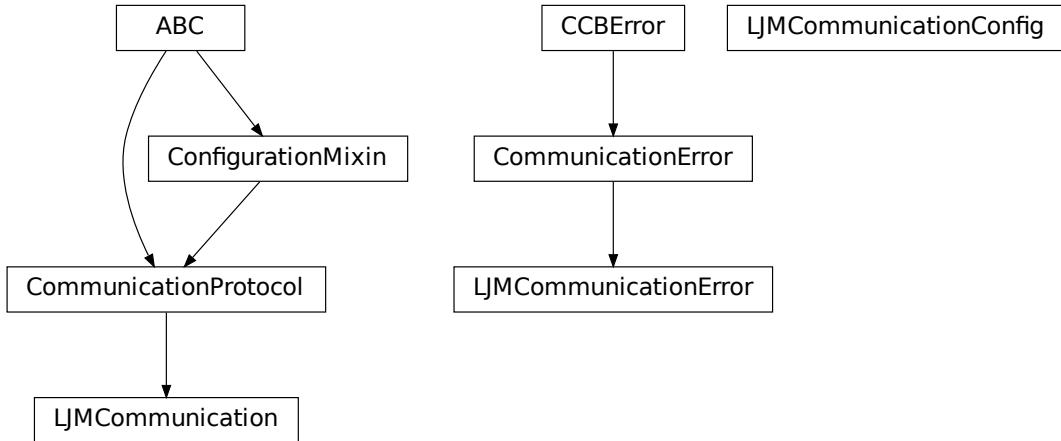

Communication protocol for LabJack using the LJM Library. Originally developed and tested for LabJack T7-PRO.

Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

class LJMCommunication(configuration)

Bases: `CommunicationProtocol`

Communication protocol implementing the LabJack LJM Library Python wrapper.

close() → None

Close the communication port.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property is_open: bool

Flag indicating if the communication port is open.

Returns

`True` if the port is open, otherwise `False`

open() → None

Open the communication port.

read_name(*names: str, return_num_type: typefromnumbers.Real] = <class 'float'>) → Real | Sequence[Real]

Read one or more input numeric values by name.

Parameters

- **names** – one or more names to read out from the LabJack
- **return_num_type** – optional numeric type specification for return values; by default `float`.

Returns

answer of the LabJack, either single number or multiple numbers in a sequence, respectively, when one or multiple names to read were given

Raises

TypeError – if read value of type not compatible with *return_num_type*

write_name(*name*: str, *value*: Real) → *None*

Write one value to a named output.

Parameters

- **name** – String or with name of LabJack IO
- **value** – is the value to write to the named IO port

write_names(*name_value_dict*: dict[str, Real]) → *None*

Write more than one value at once to named outputs.

Parameters

name_value_dict – is a dictionary with string names of LabJack IO as keys and corresponding numeric values

class LJMCommunicationConfig(*device_type*: str | DeviceType = 'ANY', *connection_type*: str | ConnectionType = 'ANY', *identifier*: str = 'ANY')

Bases: object

Configuration dataclass for [LJMCommunication](#).

class ConnectionType(*value*=<no_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: [AutoNumberNameEnum](#)

LabJack connection type.

ANY = 1

ETHERNET = 4

TCP = 3

USB = 2

WIFI = 5

class DeviceType(*value*=<no_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: [AutoNumberNameEnum](#)

LabJack device types.

Can be also looked up by ambiguous Product ID (*p_id*) or by instance name: `python LabJackDeviceType(4)` is `LabJackDeviceType('T4')`

ANY = 1

T4 = 2

T7 = 3

T7_PRO = 4

classmethod `get_by_p_id(p_id: int) → DeviceType | list[DeviceType]`

Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

Parameters

`p_id` – Product ID of a LabJack device

Returns

 Instance or list of instances of `LabJackDeviceType`

Raises

`ValueError` – when Product ID is unknown

clean_values() → `None`

Performs value checks on `device_type` and `connection_type`.

connection_type: `str | ConnectionType = 'ANY'`

Can be either string or of enum `ConnectionType`.

device_type: `str | DeviceType = 'ANY'`

Can be either string ‘ANY’, ‘T7_PRO’, ‘T7’, ‘T4’, or of enum `DeviceType`.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- `fieldname` – name of the field
- `value` – value to assign

identifier: `str = 'ANY'`

The identifier specifies information for the connection to be used. This can be an IP address, serial number, or device name. See the LabJack docs (<https://labjack.com/support/software/api/ljm/function-reference/ljmopens/identifier-parameter>) for more information.

is_configdataclass = True

classmethod keys() → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

Returns

 a list of strings containing all keys.

classmethod optional_defaults() → `dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

 a list of strings containing all optional keys.

classmethod required_keys() → `Sequence[str]`


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

 a list of strings containing all required keys.

exception LJMCommunicationErrorBases: *CommunicationError*

Errors coming from LJMCommunication.

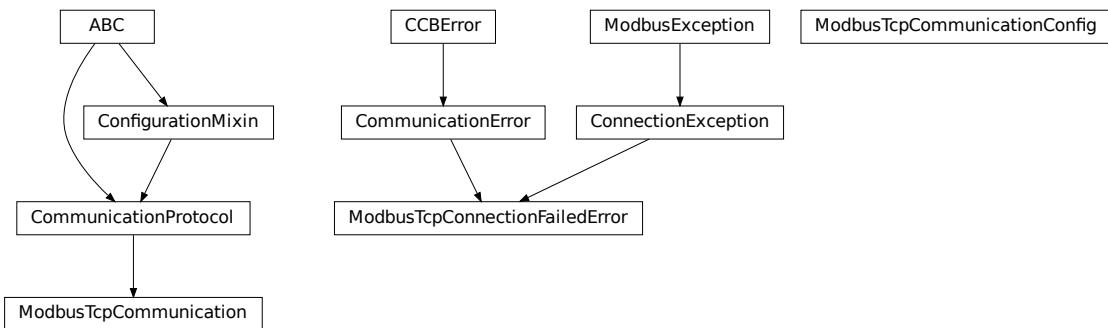hvl_ccb.comm.modbus_tcpCommunication protocol for modbus TCP ports. Makes use of the `pymodbus` library.**class ModbusTcpCommunication(configuration)**Bases: *CommunicationProtocol*

Implements the Communication Protocol for modbus TCP.

close()

Close the Modbus TCP connection.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

open() → None

Open the Modbus TCP connection.

Raises`ModbusTcpConnectionFailedError` – if the connection fails.**read_holding_registers(address: int, count: int) → list[int]**

Read specified number of register starting with given address and return the values from each register.

Parameters

- **address** – address of the first register
- **count** – count of registers to read

Returnslist of `int` values**read_input_registers(address: int, count: int) → list[int]**

Read specified number of register starting with given address and return the values from each register in a list.

Parameters

- **address** – address of the first register
- **count** – count of registers to read

Returns

list of *int* values

write_registers(*address*: *int*, *values*: *list[int]* | *int*)

Write values from the specified address forward.

Parameters

- **address** – address of the first register
- **values** – list with all values

class ModbusTcpCommunicationConfig(*host*: *str* | *IPv4Address* | *IPv6Address*, *unit*: *int*, *port*: *int* = 502)

Bases: *object*

Configuration dataclass for *ModbusTcpCommunication*.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: *str* | *IPv4Address* | *IPv6Address*

is_configdataclass = *True*

classmethod keys() → *Sequence[str]*

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → *dict[str, object]*

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

port: *int* = 502

classmethod required_keys() → *Sequence[str]*


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
unit: int

exception ModbusTcpConnectionFailedError(string= "")
    Bases: ConnectionException, CommunicationError
    Error raised when the connection failed.
```

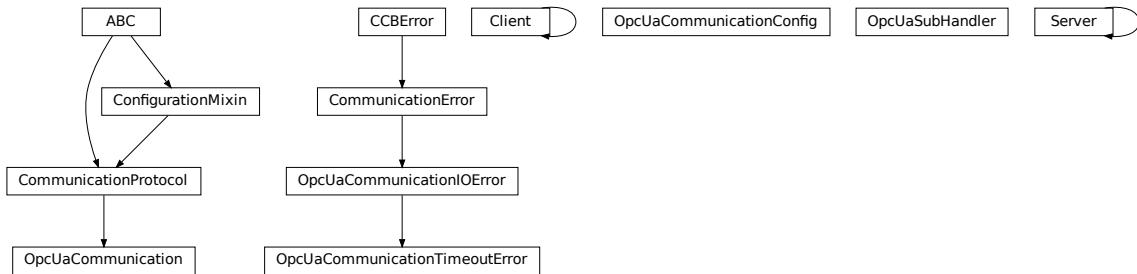hvl_ccb.comm.opc

Communication protocol implementing an OPC UA connection. This protocol is used to interface with the “Cube” PLC from Siemens.

```
class Client(url: str, timeout: int = 4)
    Bases: Client
    disconnect()
    get_objects_node()
        Get Objects node of client. Returns a Node object.
    property is_open
    send_hello(*args, **kwargs)

class OpcUaCommunication(config)
    Bases: CommunicationProtocol
    Communication protocol implementing an OPC UA connection. Makes use of the package python-opcua.
    close() → None
        Close the connection to the OPC UA server.
    static config_cls()
        Return the default configdataclass class.

        Returns
            a reference to the default configdataclass class

    init_monitored_nodes(node_id: object | Iterable, ns_index: int) → None
        Initialize monitored nodes.

        Parameters
            • node_id – one or more strings of node IDs; node IDs are always casted via str() method here, hence do not have to be strictly string objects.
```

- **ns_index** – the namespace index the nodes belong to.

Raises

OpcUaCommunicationIOError – when protocol was not opened or can't communicate with a OPC UA server

property is_open: bool

Flag indicating if the communication port is open. —DEPRECATED! DO NOT USE!!!—

Returns

True if the port is open, otherwise *False*

open() → None

Open the communication to the OPC UA server.

Raises

OpcUaCommunicationIOError – when communication port cannot be opened.

read(node_id, ns_index)

Read a value from a node with id and namespace index.

Parameters

- **node_id** – the ID of the node to read the value from
- **ns_index** – the namespace index of the node

Returns

the value of the node object.

Raises

OpcUaCommunicationIOError – when protocol was not opened or can't communicate with a OPC UA server

write(node_id, ns_index, value) → None

Write a value to a node with name **name**.

Parameters

- **node_id** – the id of the node to write the value to.
- **ns_index** – the namespace index of the node.
- **value** – the value to write.

Raises

OpcUaCommunicationIOError – when protocol was not opened or can't communicate with a OPC UA server

```
class OpcUaCommunicationConfig(host: str | ~ipaddress.IPv4Address | ~ipaddress.IPv6Address,
                                endpoint_name: str, port: int = 4840, sub_handler:
                                ~hvl_ccb.comm.opc.OpcUaSubHandler =
                                <hvl_ccb.comm.opc.OpcUaSubHandler object>, update_parameter:
                                ~asyncua.ua.uaproto.protocol_auto.CreateSubscriptionParameters =
                                CreateSubscriptionParameters(RequestedPublishingInterval=1000,
                                RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22,
                                MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0),
                                wait_timeout_retry_sec: int | float = 1, max_timeout_retry_nr: int = 5)
```

Bases: **object**

Configuration dataclass for OPC UA Communciation.

clean_values()

endpoint_name: str

Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str | IPv4Address | IPv6Address

Hostname or IP-Address of the OPC UA server.

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

max_timeout_retry_nr: int = 5

Maximal number of call re-tries on underlying OPC UA client timeout error

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

port: int = 4840

Port of the OPC UA server to connect to.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

sub_handler: OpcUaSubHandler = <hvl_ccb.comm.opc.OpcUaSubHandler object>

object to use for handling subscriptions.

update_parameter: CreateSubscriptionParameters = CreateSubscriptionParameters(RequestedPublishingInterval=1000, RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22, MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0)

Values are given as a *ua.CreateSubscriptionParameters* as these parameters are requested by the OPC server. Other values will lead to an automatic revision of the parameters and a warning in the opc-logger, cf. MR !173

```
wait_timeout_retry_sec: int | float = 1
    Wait time between re-trying calls on underlying OPC UA client timeout error

exception OpcUaCommunicationIOError
    Bases: OSError, CommunicationError
    OPC-UA communication I/O error.

exception OpcUaCommunicationTimeoutError
    Bases: OpcUaCommunicationIOError
    OPC-UA communication timeout error.


class OpcUaSubHandler
    Bases: object
    Base class for subscription handling of OPC events and data change events. Override methods from this class to add own handling capabilities.

    To receive events from server for a subscription data_change and event methods are called directly from receiving thread. Do not do expensive, slow or network operation there. Create another thread if you need to do such a thing.

    datachange_notification(node, val, data)
    event_notification(event)

class Server(shelf_file=None, tloop=None, sync_wrapper_timeout: float | None = 120)
    Bases: Server
    get_objects_node()
        Get Objects node of server. Returns a Node object.
```

hvl_ccb.comm.serial

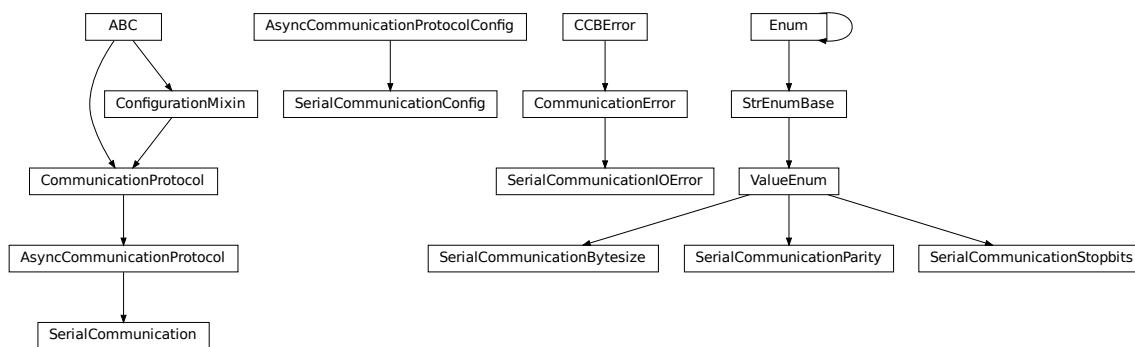

Communication protocol for serial ports. Makes use of the `pySerial` library.

```
class SerialCommunication(configuration)
    Bases: AsyncCommunicationProtocol
    Implements the Communication Protocol for serial ports.
```

close()

Close the serial connection.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property is_open: bool

Flag indicating if the serial port is open.

Returns

True if the serial port is open, otherwise *False*

open()

Open the serial connection.

Raises

SerialCommunicationIOError – when communication port cannot be opened.

read_bytes() → bytes

Read the bytes from the serial port till the terminator is found. The input buffer may hold additional lines afterwards.

This method uses *self.access_lock* to ensure thread-safety.

Returns

Bytes read from the serial port; *b”* if there was nothing to read.

Raises

SerialCommunicationIOError – when communication port is not opened

read_single_bytes(size: int = 1) → bytes

Read the specified number of bytes from the serial port. The input buffer may hold additional data afterwards.

Returns

Bytes read from the serial port; *b”* if there was nothing to read.

write_bytes(data: bytes) → int

Write bytes to the serial port.

This method uses *self.access_lock* to ensure thread-safety.

Parameters

data – data to write to the serial port

Returns

number of bytes written

Raises

SerialCommunicationIOError – when communication port is not opened

class SerialCommunicationBytesize(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)

Bases: *ValueEnum*

Serial communication bytesize.

EIGHTBITS = 8

```
FIVEBITS = 5
SEVENBITS = 7
SIXBITS = 6

class SerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: int
                                | float = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: str |
                                None = None, baudrate: int = 9600, parity: str |
                                SerialCommunicationParity = SerialCommunicationParity.NONE,
                                stopbits: int | float | SerialCommunicationStopbits =
                                SerialCommunicationStopbits.ONE, bytesize: int |
                                SerialCommunicationBytesize =
                                SerialCommunicationBytesize.EIGHTBITS, timeout: int | float = 2)
```

Bases: [AsyncCommunicationProtocolConfig](#)

Configuration dataclass for [SerialCommunication](#).

Bytesize

alias of [SerialCommunicationBytesize](#)

Parity

alias of [SerialCommunicationParity](#)

Stopbits

alias of [SerialCommunicationStopbits](#)

baudrate: int = 9600

Baudrate of the serial port

bytesize: int | SerialCommunicationBytesize = 8

Size of a byte, 5 to 8

clean_values()

create_serial_port() → Serial

Create a serial port instance according to specification in this configuration

Returns

Closed serial port instance

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

```
classmethod optional_defaults() → dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

```
parity: str | SerialCommunicationParity = 'N'
```

Parity to be used for the connection.

```
port: str | None = None
```

Port is a string referring to a COM-port (e.g. 'COM3') or a URL. The full list of capabilities is found [on the pyserial documentation](#).

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
stopbits: int | float | SerialCommunicationStopbits = 1
```

Stopbits setting, can be 1, 1.5 or 2.

```
terminator_str() → str
```

```
timeout: int | float = 2
```

Timeout in seconds for the serial port

```
exception SerialCommunicationIOError
```

Bases: OSError, [CommunicationError](#)

Serial communication related I/O errors.

```
class SerialCommunicationParity(value=<no_arg>, names=None, module=None, qualname=None,  
                                 type=None, start=1, boundary=None)
```

Bases: [ValueEnum](#)

Serial communication parity.

```
EVEN = 'E'
```

```
MARK = 'M'
```

```
NAMES = {'E': 'Even', 'M': 'Mark', 'N': 'None', 'O': 'Odd', 'S': 'Space'}
```

```
NONE = 'N'
```

```
ODD = 'O'
```

```
SPACE = 'S'
```

```
class SerialCommunicationStopbits(value=<no_arg>, names=None, module=None, qualname=None,  
                                   type=None, start=1, boundary=None)
```

Bases: [ValueEnum](#)


Serial communication stopbits.

```
ONE = 1
```

```
ONE_POINT_FIVE = 1.5
```

```
TWO = 2
```

hvl_ccb.comm.tcp

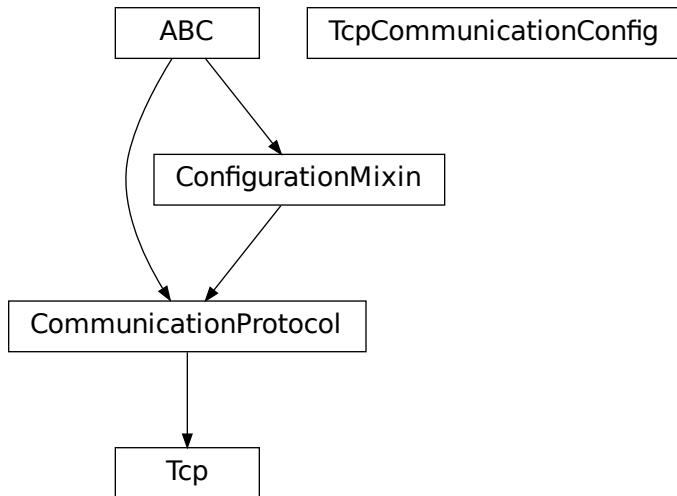

Makes use of the socket library.

TCP communication protocol.

class `Tcp(configuration)`

Bases: `CommunicationProtocol`

Tcp Communication Protocol.

close() → `None`

Close TCP connection.

static config_cls() → type[`TcpCommunicationConfig`]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

open() → `None`

Open TCP connection.

read() → str

TCP read function :return: information read from TCP buffer formatted as string

write(command: str = "") → `None`

TCP write function :param command: command string to be sent :return: none

```
class TcpCommunicationConfig(host: str | IPv4Address | IPv6Address, port: int = 54321, bufsize: int = 1024)
Bases: object
Configuration dataclass for TcpCommunication.

bufsize: int = 1024

clean_values()

force_value(fieldname, value)
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define post_force_value method with same signature as this method to do extra processing
after value has been forced on fieldname.

Parameters

- fieldname – name of the field
- value – value to assign


host: str | IPv4Address | IPv6Address

is_configdataclass = True

classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns
a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
on instantiation.


Returns
a list of strings containing all optional keys.

port: int = 54321

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
instantiation.

Returns
a list of strings containing all required keys.
```

hvl_ccb.comm.telnet

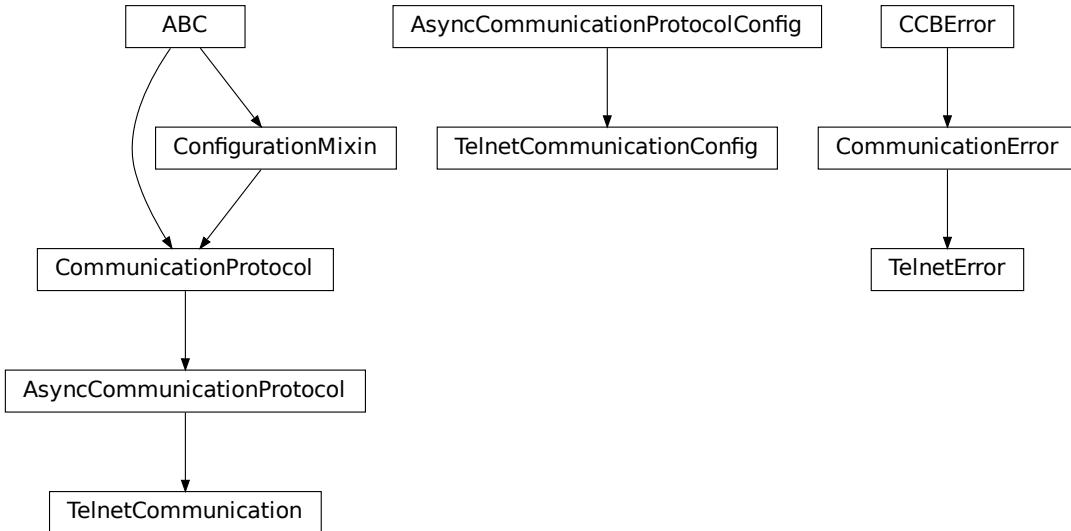

Communication protocol for telnet. Makes use of the `telnetlib` library.

class TelnetCommunication(configuration)

Bases: `AsyncCommunicationProtocol`

Implements the Communication Protocol for telnet.

close()

Close the telnet connection unless it is not closed.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property is_open: bool

Is the connection open?

Returns

True for an open connection

open()

Open the telnet connection unless it is not yet opened.

read_bytes() → bytes

Read data as *bytes* from the telnet connection.

Returns

data from telnet connection

Raises

`TelnetError` – when connection is not open, raises an Error during the communication

write_bytes(*data*: bytes)

Write the data as *bytes* to the telnet connection.

Parameters

- **data** – Data to be sent.

Raises

- **TelnetError** – when connection is not open, raises an Error during the communication

```
class TelnetCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: int
                                | float = 0.5, default_n_attempts_read_text_nonempty: int = 10, host: str |
                                IPv4Address | IPv6Address | None = None, port: int = 0, timeout: int |
                                float = 0.2)
```

Bases: *AsyncCommunicationProtocolConfig*

Configuration dataclass for *TelnetCommunication*.

clean_values()**create_telnet()** → Telnet | *None*

Create a telnet client :return: Opened Telnet object or None if connection is not possible

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str | IPv4Address | IPv6Address | *None* = *None*

Host to connect to can be localhost or

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

port: int = 0

Port at which the host is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
timeout: int | float = 0.2
```


Timeout for reading a line

exception TelnetError

Bases: `OSError`, `CommunicationError`

Telnet communication related errors.

hvl_ccb.comm.visa

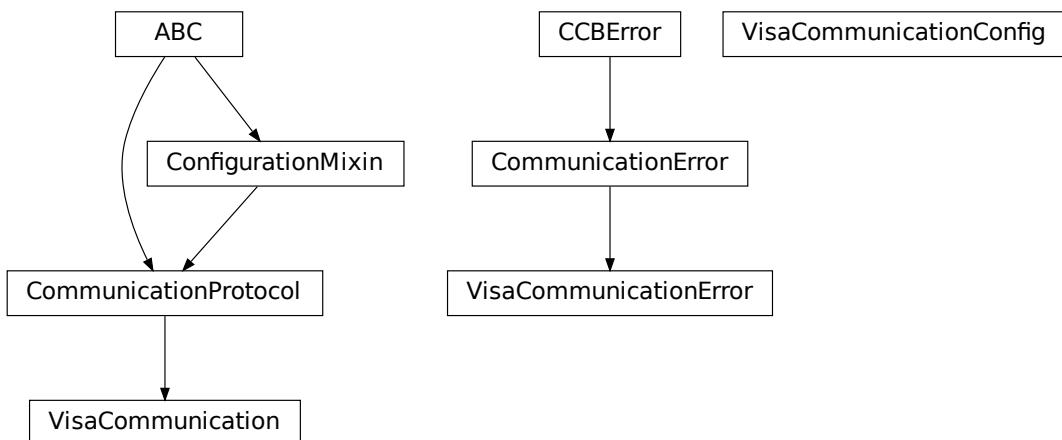

Communication protocol for VISA. Makes use of the pyvisa library. The backend can be NI-Visa or pyvisa-py.

Information on how to install a VISA backend can be found here: https://pyvisa.readthedocs.io/en/master/getting_nivisa.html

So far only TCPIP SOCKET and TCPIP INSTR interfaces are supported.

class VisaCommunication(configuration)

Bases: `CommunicationProtocol`

Implements the Communication Protocol for VISA / SCPI.

MULTI_COMMANDS_MAX = 5

The maximum of commands that can be sent in one round is 5 according to the VISA standard.

MULTI_COMMANDS_SEPARATOR = ';'

The character to separate two commands is ; according to the VISA standard.

WAIT_AFTER_WRITE = 0.08

Small pause in seconds to wait after write operations, allowing devices to really do what we tell them before continuing with further tasks.

close() → None

Close the VISA connection and invalidates the handle.

static config_cls() → type[*VisaCommunicationConfig*]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

open() → *None*

Open the VISA connection and create the resource.

query(*commands: str) → str | tuple[str, ...]

A combination of write(message) and read.

Parameters

commands – list of commands

Returns

list of values

Raises

VisaCommunicationError – when connection was not started, or when trying to issue too many commands at once.

spoll() → int

Execute serial poll on the device. Reads the status byte register STB. This is a fast function that can be executed periodically in a polling fashion.

Returns

integer representation of the status byte

Raises

VisaCommunicationError – when connection was not started

write(*commands: str) → *None*

Write commands. No answer is read or expected.

Parameters

commands – one or more commands to send

Raises

VisaCommunicationError – when connection was not started

class VisaCommunicationConfig(host: str | IPv4Address | IPv6Address, interface_type: str | InterfaceType, board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = '')

Bases: *object*

VisaCommunication configuration dataclass.

class InterfaceType(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)

Bases: *AutoNumberNameEnum*

Supported VISA Interface types.

TCPIP_INSTR = 2

VXI-11 protocol

TCPIP_SOCKET = 1

VISA-Raw protocol

address(*host*: str, *port*: int | None = None, *board*: int | None = None) → str

Address string specific to the VISA interface type.

Parameters

- **host** – host IP address
- **port** – optional TCP port
- **board** – optional board number

Returns

address string

property address: str

Address string depending on the VISA protocol's configuration.

Returns

address string corresponding to current configuration

board: int = 0

Board number is typically 0 and comes from old bus systems.

chunk_size: int = 204800

Chunk size is the allocated memory for read operations. The standard is 20kB, and is increased per default here to 200kB. It is specified in bytes.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str | IPv4Address | IPv6Address

interface_type: str | InterfaceType

Interface type of the VISA connection, being one of *InterfaceType*.

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

open_timeout: int = 1000

Timeout for opening the connection, in milli seconds.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

```

port: int = 5025
    TCP port, standard is 5025.

read_termination: str = '\n'
    Read termination character.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns
    a list of strings containing all required keys.

timeout: int = 5000
    Timeout for commands in milli seconds.

visa_backend: str = ''
    Specifies the path to the library to be used with PyVISA as a backend. Defaults to None, which is NI-VISA (if installed), or pyvisa-py (if NI-VISA is not found). To force the use of pyvisa-py, specify '@py' here.

write_termination: str = '\n'
    Write termination character.

exception VisaCommunicationError
    Bases: OSError, CommunicationError
    Base class for VisaCommunication errors.

```

Module contents

Communication protocols subpackage.


hvl_ccb.dev

Subpackages

hvl_ccb.dev.crylas

Submodules

hvl_ccb.dev.crylas.crylas

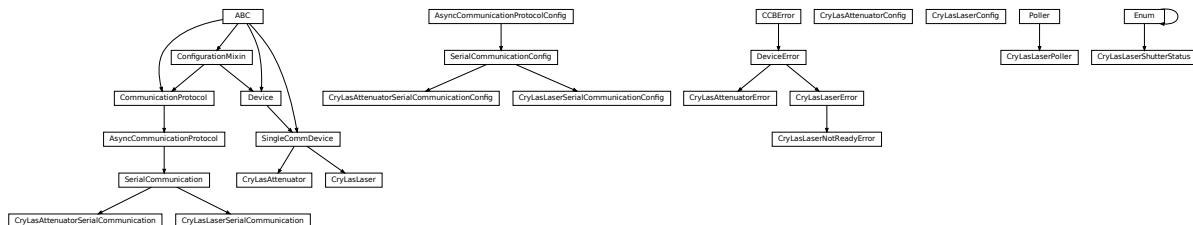

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: https://www.crylas.de/products/pulsed_laser.html

class CryLasAttenuator(*com, dev_config=None*)

Bases: *SingleCommDevice*

Device class for the CryLas laser attenuator.

property attenuation: int | float

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

set_attenuation(percent: int | float) → None

Set the percentage of attenuated light (inverse of set_transmission). :param percent: percentage of attenuation, number between 0 and 100 :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

set_init_attenuation()

Sets the attenuation to its configured initial/default value

Raises

SerialCommunicationIOError – when communication port is not opened

set_transmission(percent: int | float) → None

Set the percentage of transmitted light (inverse of set_attenuation). :param percent: percentage of transmitted light :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

start() → None

Open the com, apply the config value ‘init_attenuation’

Raises

SerialCommunicationIOError – when communication port cannot be opened

property transmission: int | float

class CryLasAttenuatorConfig(*init_attenuation: int | float = 0, response_sleep_time: int | float = 1*)

Bases: *object*

Device configuration dataclass for CryLas attenuator.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

init_attenuation: int | float = 0**is_configdataclass** = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

response_sleep_time: int | float = 1**exception CryLasAttenuatorError**

Bases: *DeviceError*

General error with the CryLas Attenuator.

class CryLasAttenuatorSerialCommunication(*configuration*)

Bases: *SerialCommunication*

Specific communication protocol implementation for the CryLas attenuator. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class CryLasAttenuatorSerialCommunicationConfig(terminator: bytes = b'', encoding: str = 'utf-8',
                                                encoding_error_handling: str = 'strict',
                                                wait_sec_read_text_nonempty: Union[int, float] =
                                                0.5, default_n_attempts_read_text_nonempty: int =
                                                10, port: Optional[str] = None, baudrate: int = 9600,
                                                parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 3)
```

Bases: *SerialCommunicationConfig*

baudrate: int = 9600

Baudrate for CryLas attenuator is 9600 baud

bytesize: int | SerialCommunicationBytesize = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

CryLas attenuator does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```

stopbits: int | SerialCommunicationStopbits = 1
    CryLas attenuator uses one stop bit

terminator: bytes = b''
    No terminator

timeout: int | float = 3
    use 3 seconds timeout as default

class CryLasLaser(com, dev_config=None)
    Bases: SingleCommDevice

    CryLas laser controller device class.

class AnswersShutter(value=<no_arg>, names=None, module=None, qualname=None, type=None,
start=1, boundary=None)
    Bases: Enum

    Standard answers of the CryLas laser controller to 'Shutter' command passed via com.

    CLOSED = 'Shutter inaktiv'

    OPENED = 'Shutter aktiv'

class AnswersStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None,
start=1, boundary=None)
    Bases: Enum

    Standard answers of the CryLas laser controller to 'STATUS' command passed via com.

    ACTIVE = 'STATUS: Laser active'

    HEAD = 'STATUS: Head ok'

    INACTIVE = 'STATUS: Laser inactive'

    READY = 'STATUS: System ready'

    TEC1 = 'STATUS: TEC1 Regulation ok'

    TEC2 = 'STATUS: TEC2 Regulation ok'

class LaserStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
boundary=None)
    Bases: Enum

    Status of the CryLas laser

    READY_ACTIVE = 2

    READY_INACTIVE = 1

    UNREADY_INACTIVE = 0

    property is_inactive

    property is_ready

```

```
class RepetitionRates(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                      start=1, boundary=None)
```

Bases: IntEnum

Repetition rates for the internal software trigger in Hz

HARDWARE = 0

SOFTWARE_INTERNAL_SIXTY = 60

SOFTWARE_INTERNAL_TEN = 10

SOFTWARE_INTERNAL_TWENTY = 20

ShutterStatus

alias of [CryLasLaserShutterStatus](#)

close_shutter() → *None*

Close the laser shutter.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

get_pulse_energy_and_rate() → tuple[int, int]

Use the debug mode, return the measured pulse energy and rate.

Returns

(energy in micro joule, rate in Hz)

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if the device does not answer the query

laser_off() → *None*

Turn the laser off.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

laser_on() → *None*

Turn the laser on.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserNotReadyError** – if the laser is not ready to be turned on
- **CryLasLaserError** – if success is not confirmed by the device

open_shutter() → *None*

Open the laser shutter.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

set_init_shutter_status() → *None*

Open or close the shutter, to match the configured shutter_status.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

set_pulse_energy(energy: int) → *None*

Sets the energy of pulses (works only with external hardware trigger). Proceed with small energy steps, or the regulation may fail.

Parameters

energy – energy in micro joule

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if the device does not confirm success

set_repetition_rate(rate: int | RepetitionRates) → *None*

Sets the repetition rate of the internal software trigger.

Parameters

rate – frequency (Hz) as an integer

Raises

- **ValueError** – if rate is not an accepted value in RepetitionRates Enum
- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

start() → *None*

Opens the communication protocol and configures the device.

Raises

SerialCommunicationIOError – when communication port cannot be opened

stop() → *None*

Stops the device and closes the communication protocol.

Raises

- **SerialCommunicationIOError** – if com port is closed unexpectedly
- **CryLasLaserError** – if laser_off() or close_shutter() fail

```
property target_pulse_energy

update_laser_status() → None
    Update the laser status to LaserStatus.NOT_READY or LaserStatus.INACTIVE or LaserStatus.ACTIVE.
    Note: laser never explicitly says that it is not ready (LaserStatus.NOT_READY) in response to 'STATUS' command. It only says that it is ready (heated-up and implicitly inactive/off) or active (on). If it's not either of these then the answer is Answers.HEAD. Moreover, the only time the laser explicitly says that its status is inactive (Answers.INACTIVE) is after issuing a 'LASER OFF' command.

    Raises
        SerialCommunicationIOError – when communication port is not opened

update_repetition_rate() → None
    Query the laser repetition rate.

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • CryLasLaserError – if success is not confirmed by the device

update_shutter_status() → None
    Update the shutter status (OPENED or CLOSED)

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • CryLasLaserError – if success is not confirmed by the device

update_target_pulse_energy() → None
    Query the laser pulse energy.

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • CryLasLaserError – if success is not confirmed by the device

wait_until_ready() → None
    Block execution until the laser is ready

    Raises
        CryLasLaserError – if the polling thread stops before the laser is ready

class CryLasLaserConfig(calibration_factor: int | float = 4.35, polling_period: int | float = 12, polling_timeout: int | float = 300, auto_laser_on: bool = True, init_shutter_status: int | CryLasLaserShutterStatus = CryLasLaserShutterStatus.CLOSED)
    Bases: object
    Device configuration dataclass for the CryLas laser controller.

    ShutterStatus
        alias of CryLasLaserShutterStatus
        auto_laser_on: bool = True
        calibration_factor: int | float = 4.35
        clean_values()
```

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

init_shutter_status: int | *CryLasLaserShutterStatus* = 0**is_configdataclass** = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

polling_period: int | float = 12**polling_timeout:** int | float = 300**classmethod required_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

exception CryLasLaserError

Bases: *DeviceError*

General error with the CryLas Laser.

exception CryLasLaserNotReadyError

Bases: *CryLasLaserError*

Error when trying to turn on the CryLas Laser before it is ready.

class CryLasLaserPoller(*spoll_handler*: Callable, *check_handler*: Callable, *check_laser_status_handler*: Callable, *polling_delay_sec*: int | float = 0, *polling_interval_sec*: int | float = 1, *polling_timeout_sec*: int | float | None = None)

Bases: *Poller*

Poller class for polling the laser status until the laser is ready.

Raises

- **CryLasLaserError** – if the timeout is reached before the laser is ready
- **SerialCommunicationIOError** – when communication port is closed.

class CryLasLaserSerialCommunication(*configuration*)

Bases: *SerialCommunication*

Specific communication protocol implementation for the CryLas laser controller. Already predefines device-specific protocol parameters in config.

READ_TEXT_SKIP_PREFIXES = ('>', 'MODE:')

Prefixes of lines that are skipped when read from the serial port.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

query(*cmd*: str, *prefix*: str, *post_cmd*: str | None = None) → str

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns the line in question.

Parameters

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer
- **post_cmd** – optional additional command to send after the query

Returns

line in question as a string

Raises

SerialCommunicationIOError – when communication port is not opened

query_all(*cmd*: str, *prefix*: str)

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns a list of successive lines starting with prefix.

Parameters

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer

Returns

line in question as a string

Raises

SerialCommunicationIOError – when communication port is not opened

read() → str

Read first line of text from the serial port that does not start with any of *self.READ_TEXT_SKIP_PREFIXES*.

Returns

String read from the serial port; '' if there was nothing to read.

Raises

SerialCommunicationIOError – when communication port is not opened

```
class CryLasLaserSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Optional[str] = None, baudrate: int = 19200, parity:
                                             Union[str,
                                               hvl_ccb.comm.serial.SerialCommunicationParity] =
                                               <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                               Union[int,
                                                 hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                 <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                 Union[int,
                                                   hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                   <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                                   Union[int, float] = 10)
```

Bases: *SerialCommunicationConfig*

baudrate: int = 19200

Baudrate for CryLas laser is 19200 baud

bytesize: int | SerialCommunicationBytesize = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

CryLas laser does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
stopbits: int | SerialCommunicationStopbits = 1
    CryLas laser uses one stop bit
terminator: bytes = b'\n'
    The terminator is LF
timeout: int | float = 10
    use 10 seconds timeout as default (a long timeout is needed!)
class CryLasLaserShutterStatus(value=<no_arg>, names=None, module=None, qualname=None,
                                 type=None, start=1, boundary=None)
Bases: Enum
Status of the CryLas laser shutter
CLOSED = 0
OPENED = 1
```

Module contents

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: https://www.crylas.de/products/pulsed_laser.html


hvl_ccb.dev.cube

Submodules

hvl_ccb.dev.cube.alarms

Alarms

Alarms of the different “Cubes”.

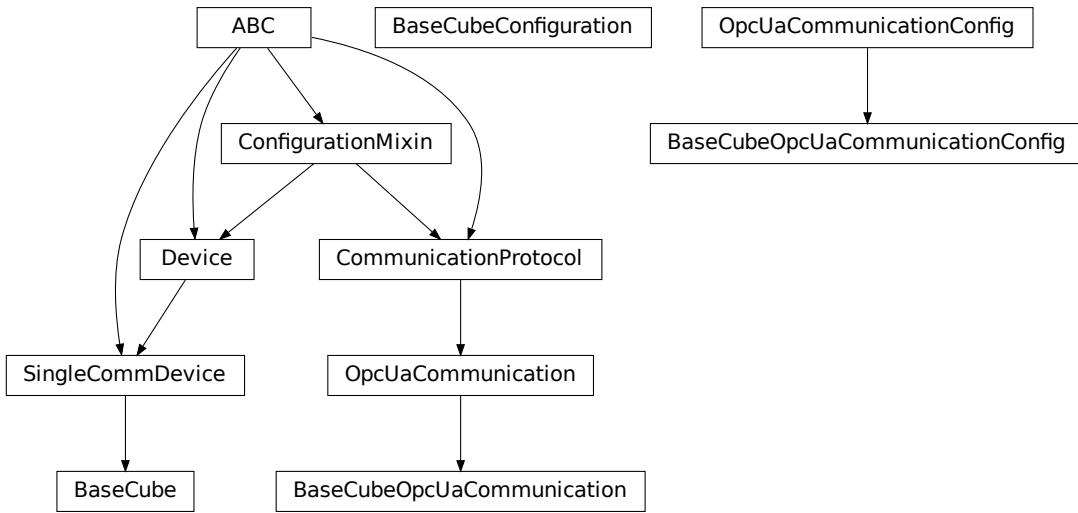hvl_ccb.dev(cube).base

Classes for the BaseCube device.

```

class BaseCube(com, dev_config=None)
  Bases: SingleCommDevice
  Base class for Cube variants.

  OPC_MAX_YEAR = 2089
  OPC_MIN_YEAR = 1990

  active_alarms(human_readable: bool = True) → list[int | str]
    Displays all active alarms / messages.

    Parameters
      human_readable – True for human readable message, False for corresponding integer

    Returns
      list with active alarms

  property breakdown_detection_active: bool
    Get the state of the breakdown detection functionality. Returns True if it is enabled, False otherwise.

    Returns
      state of the breakdown detection functionality

  breakdown_detection_reset() → None
    Reset the breakdown detection circuitry so that it is ready to detect breakdowns again.

  property breakdown_detection_triggered: bool
    See if breakdown detection unit has been triggered. Returns True if it is triggered, False otherwise.

    Returns
      trigger status of the breakdown detection unit
  
```

property cee16_socket

Read the on-state of the IEC CEE16 three-phase power socket.

Returns

the on-state of the CEE16 power socket

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

classmethod datetime_to_opc(*time_dt*: datetime) → list[int]

Converts python datetime format into opc format (list of 8 integers) as defined in the following link: <https://support.industry.siemens.com/cs/mdm/109798671?c=133950752267&lc=de-WW> Each byte corresponds to one list entry. [yy, MM, dd, hh, mm, ss, milliseconds, weekday] Milliseconds and Weekday are not used, as this precision / information is not needed. The conversion of the numbers is special. Each decimal number is treated as it would be a hex-number and then converted back to decimal. This is tested with the used PLC in the BaseCube. yy: 0 to 99 (0 -> 2000, 89 -> 2089, 90 -> 1990, 99 -> 1999) MM: 1 to 12 dd: 1 to 31 hh: 0 to 23 mm: 0 to 59 ss: 0 to 59

Parameters

time_dt – time to be converted

Returns

time in opc list format

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

display_message_board() → None

Display 15 newest messages

display_status_board() → None

Display status board.

door_1_status

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

door_2_status

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

door_3_status

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

earthing_rod_1_status

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

earthing_rod_2_status

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

earthing_rod_3_status

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

property operate: bool | None

Indicates if ‘operate’ is activated. ‘operate’ means locket safety circuit, red lamps, high voltage on and locked safety switches.

Returns

True if operate is activated (RED_OPERATE), *False* if ready is deactivated (RED_READY),
None otherwise

quit_error() → None

Quits errors that are active on the Cube.

read(node_id: str)

Local wrapper for the OPC UA communication protocol read method.

Parameters

node_id – the id of the node to read.

Returns

the value of the variable

property ready: bool | None

Indicates if ‘ready’ is activated. ‘ready’ means locket safety circuit, red lamps, but high voltage still off.

Returns

True if ready is activated (RED_READY), *False* if ready is deactivated (GREEN_READY),
None otherwise

set_message_board(msgs: list[str], display_board: bool = True) → None

Fills messages into message board that display that 15 newest messages with a timestamp.

Parameters

- **msgs** – list of strings
- **display_board** – display 15 newest messages if *True* (default)

Raises

ValueError – if there are too many messages or the positions indices are invalid.

set_status_board(msgs: list[str], pos: list[int] | None = None, clear_board: bool = True, display_board: bool = True) → None

Sets and displays a status board. The messages and the position of the message can be defined.

Parameters

- **msgs** – list of strings
- **pos** – list of integers [0...14]
- **clear_board** – clear unspecified lines if *True* (default), keep otherwise
- **display_board** – display new status board if *True* (default)

Raises

ValueError – if there are too many messages or the positions indices are invalid.

start() → None

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

property status: SafetyStatus

Get the safety circuit status of the Cube. This methods is for the user.

Returns

the safety status of the Cube's state machine.

stop() → None

Stop the Cube device. Deactivates the remote control and closes the communication protocol.

Raises

CubeStopError – when the cube is not in the correct status to stop the operation

t13_socket_1

Set and get the state of a SEV T13 power socket.

t13_socket_2

Set and get the state of a SEV T13 power socket.

t13_socket_3

Set and get the state of a SEV T13 power socket.

write(node_id, value) → None

Local wrapper for the OPC UA communication protocol write method.

Parameters

- **node_id** – the id of the node to write
- **value** – the value to write to the variable

class BaseCubeConfiguration(namespace_index: int = 3, polling_delay_sec: int | float = 5.0, polling_interval_sec: int | float = 1.0, timeout_status_change: int | float = 6, timeout_interval: int | float = 0.1, noise_level_measurement_channel_1: int | float = 100, noise_level_measurement_channel_2: int | float = 100, noise_level_measurement_channel_3: int | float = 100, noise_level_measurement_channel_4: int | float = 100)

Bases: object

Configuration dataclass for the BaseCube devices.

clean_values()

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

```

namespace_index: int = 3
    Namespace of the OPC variables, typically this is 3 (coming from Siemens)

noise_level_measurement_channel_1: int | float = 100
noise_level_measurement_channel_2: int | float = 100
noise_level_measurement_channel_3: int | float = 100
noise_level_measurement_channel_4: int | float = 100

classmethod optional_defaults() → dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns
        a list of strings containing all optional keys.

polling_delay_sec: int | float = 5.0
polling_interval_sec: int | float = 1.0

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns
        a list of strings containing all required keys.

timeout_interval: int | float = 0.1
timeout_status_change: int | float = 6

class BaseCubeOpcUaCommunication(config)
    Bases: OpcUaCommunication
    Communication protocol specification for BaseCube devices.

static config_cls()
    Return the default configdataclass class.

    Returns
        a reference to the default configdataclass class

class BaseCubeOpcUaCommunicationConfig(host: str | ~ipaddress.IPv4Address | ~ipaddress.IPv6Address,
    endpoint_name: ~hvl_ccb.dev(cube.constants._CubeOpcEndpoint
    = _CubeOpcEndpoint.BASE_CUBE, port: int = 4840,
    sub_handler: ~hvl_ccb.comm.opc.OpcUaSubHandler =
    <hvl_ccb.dev(cube.base._BaseCubeSubscriptionHandler object>,
update_parameter:
~asyncua.ua.uaproto.protocol_auto.CreateSubscriptionParameters =
CreateSubscriptionParameters(RequestedPublishingInterval=1000,
RequestedLifetimeCount=300,
RequestedMaxKeepAliveCount=22,
MaxNotificationsPerPublish=10000, PublishingEnabled=True,
Priority=0), wait_timeout_retry_sec: int | float = 1,
max_timeout_retry_nr: int = 5)

```

Bases: *OpcUaCommunicationConfig*

Communication protocol configuration for OPC UA, specifications for the BaseCube devices.

endpoint_name: *_CubeOpcEndpoint* = 'BaseCube'

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

sub_handler: *OpcUaSubHandler* = <hvl_ccb.dev.cube.base._BaseCubeSubscriptionHandler object>

Subscription handler for data change events

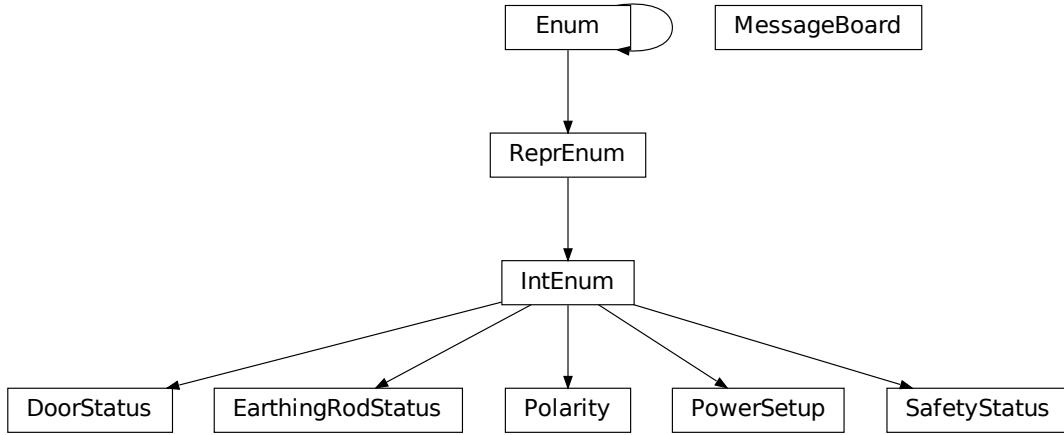hvl_ccb.dev(cube.constants)

Constants, variable names for the BaseCube OPC-connected devices.

```
class DoorStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: IntEnum

Possible status values for doors.

CLOSED = 2

Door is closed, but not locked.

ERROR = 4

Door has an error or was opened in locked state (either with emergency stop or from the inside).

INACTIVE = 0

not enabled in BaseCube HMI setup, this door is not supervised.

LOCKED = 3

Door is closed and locked (safe state).

OPEN = 1

Door is open.

```
class EarthingRodStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: IntEnum

Possible status values for earthing rods.

EXPERIMENT_BLOCKED = 0

earthing rod is somewhere in the experiment and blocks the start of the experiment

EXPERIMENT_READY = 1

earthing rod is hanging next to the door, experiment is ready to operate

```
class MessageBoard(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                    boundary=None)
```

Bases: _LineEnumBase

Variable NodeID strings for message board lines.

```
LINE_1 = '"DB_OPC_Connection"."Is_status_Line_1"'
LINE_10 = '"DB_OPC_Connection"."Is_status_Line_10"'
LINE_11 = '"DB_OPC_Connection"."Is_status_Line_11"'
LINE_12 = '"DB_OPC_Connection"."Is_status_Line_12"'
LINE_13 = '"DB_OPC_Connection"."Is_status_Line_13"'
LINE_14 = '"DB_OPC_Connection"."Is_status_Line_14"'
LINE_15 = '"DB_OPC_Connection"."Is_status_Line_15"'
LINE_2 = '"DB_OPC_Connection"."Is_status_Line_2"'
LINE_3 = '"DB_OPC_Connection"."Is_status_Line_3"'
LINE_4 = '"DB_OPC_Connection"."Is_status_Line_4"'
LINE_5 = '"DB_OPC_Connection"."Is_status_Line_5"'
LINE_6 = '"DB_OPC_Connection"."Is_status_Line_6"'
LINE_7 = '"DB_OPC_Connection"."Is_status_Line_7"'
LINE_8 = '"DB_OPC_Connection"."Is_status_Line_8"'
LINE_9 = '"DB_OPC_Connection"."Is_status_Line_9"
```

```
class Polarity(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
```

Bases: IntEnum

An enumeration.

```
NEGATIVE = 0
```

```
POSITIVE = 1
```

```
class PowerSetup(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                     boundary=None)
```

Bases: IntEnum

Possible power setups corresponding to the value of variable Power.setup. The values for slope_min are experimentally defined, below these values the slope is more like a staircase

The name of the first argument needs to be ‘value’, otherwise the IntEnum is not working correctly.

```
AC_100KV = 3
```

```
AC_150KV = 4
```

```
AC_200KV = 5
```

```
AC_50KV = 2
DC_140KV = 7
DC_280KV = 8
EXTERNAL_SOURCE = 1
IMPULSE_140KV = 9
NO_SOURCE = 0
POWER_INVERTER_220V = 6
```

STOP_SAFETY_STATUSES: tuple[*SafetyStatus*, ...] = (*SafetyStatus.GREEN_NOT_READY*,
SafetyStatus.GREEN_READY)

BaseCube's safety statuses required to close the connection to the device.

class SafetyStatus(*value=<no_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*,
boundary=None)

Bases: IntEnum

Safety status values that are possible states returned from *hvl_ccb.dev.cube.base.BaseCube.status()*.
These values correspond to the states of the BaseCube's safety circuit statemachine.

ERROR = 6

GREEN_NOT_READY = 1

GREEN_READY = 2

INITIALIZING = 0


QUICK_STOP = 5

RED_OPERATE = 4

RED_READY = 3

hvl_ccb.dev.cube.earthing_stick

EarthingStick of the different “Cubes”.

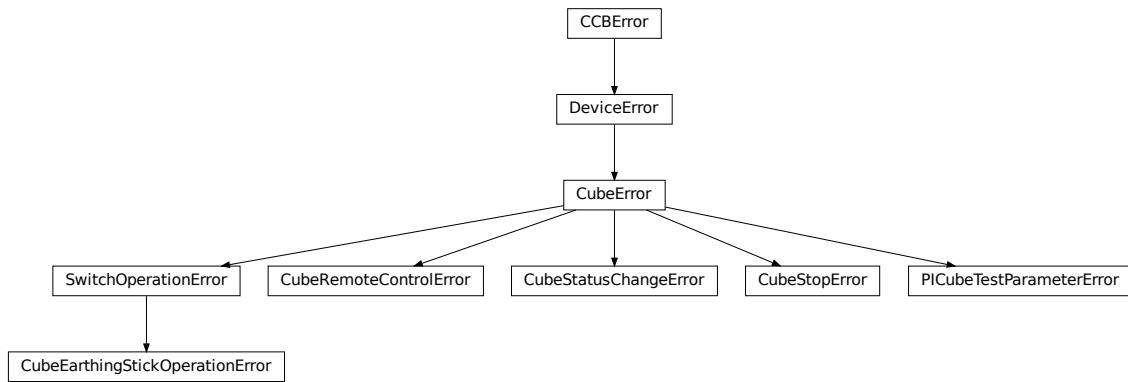hvl_ccb.dev(cube.errors

Errors of the different “Cubes”.

exception CubeEarthingStickOperationError

Bases: *SwitchOperationError*

exception CubeError

Bases: *DeviceError*

exception CubeRemoteControlError

Bases: *CubeError*

exception CubeStatusChangeError

Bases: *CubeError*

exception CubeStopError


Bases: *CubeError*

exception PICubeTestParameterError

Bases: *CubeError*

exception SwitchOperationError

Bases: *CubeError*

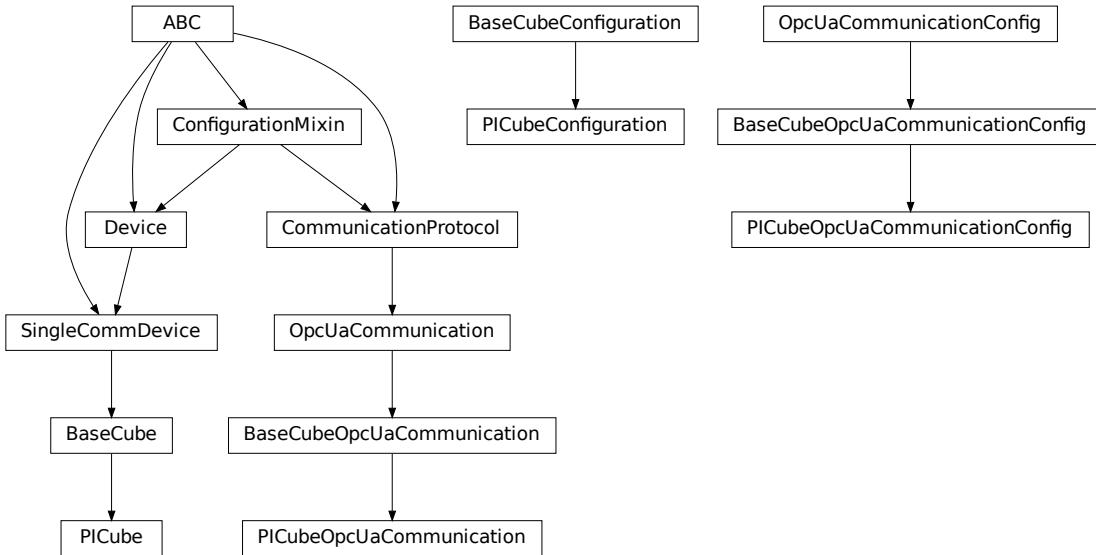hvl_ccb.dev.cube.picube

A PICube is a BaseCube with build in Power Inverter

class PICube(com, dev_config=None)

Bases: *BaseCube*

Variant of the BaseCube with build in Power Inverter

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property current_primary: float

Read the current primary current at the output of the frequency converter (before transformer).

Returns

primary current in A

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

property frequency: float

Read the electrical frequency of the current PICube setup.

Returns

the frequency in Hz

property operate: bool | None

Indicates if ‘operate’ is activated. ‘operate’ means locket safety circuit, red lamps, high voltage on and locked safety switches.

Returns

True if operate is activated (RED_OPERATE), *False* if ready is deactivated (RED_READY),
None otherwise

property polarity: Polarity | None

Polarity of a DC setup. :return: if a DC setup is programmed the polarity is returned, else None.

property power_setup: PowerSetup

Return the power setup selected in the PICube's settings.

Returns

the power setup

property voltage_actual: float

Reads the actual measured voltage and returns the value in V.

Returns

the actual voltage of the setup in V.

property voltage_max: float

Reads the maximum voltage of the setup and returns in V.

Returns

the maximum voltage of the setup in V.

property voltage_primary: float

Read the current primary voltage at the output of the frequency converter (before transformer).

Returns

primary voltage in V

```
class PICubeConfiguration(namespace_index: int = 3, polling_delay_sec: int | float = 5.0,
                         polling_interval_sec: int | float = 1.0, timeout_status_change: int | float = 6,
                         timeout_interval: int | float = 0.1, noise_level_measurement_channel_1: int | float
                         = 100, noise_level_measurement_channel_2: int | float = 100,
                         noise_level_measurement_channel_3: int | float = 100,
                         noise_level_measurement_channel_4: int | float = 100, timeout_test_parameters:
                         'Number' = 2.0)
```

Bases: *BaseCubeConfiguration*

clean_values()**force_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

```
classmethod optional_defaults() → dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
timeout_test_parameters: int | float = 2.0
```

```
class PICubeOpcUaCommunication(config)
```

Bases: *BaseCubeOpcUaCommunication*

```
static config_cls()
```

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class PICubeOpcUaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],
endpoint_name: '_CubeOpcEndpoint' =
<_CubeOpcEndpoint.PI_CUBE: 'PICube'>, port: int = 4840,
sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
<hvl_ccb.dev.cube.base._BaseCubeSubscriptionHandler object at
0x7f68981e4c10>, update_parameter:
asyncua.ua.uaproto_protocol_auto.CreateSubscriptionParameters =
CreateSubscriptionParameters(RequestedPublishingInterval=1000,
RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22,
MaxNotificationsPerPublish=10000, PublishingEnabled=True,
Priority=0), wait_timeout_retry_sec: Union[int, float] = 1,
max_timeout_retry_nr: int = 5)
```

Bases: *BaseCubeOpcUaCommunicationConfig*

```
endpoint_name: '_CubeOpcEndpoint' = 'PICube'
```

Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns

a list of strings containing all required keys.

hvl_ccb.dev(cube).support

Supports of the different “Cubes”.

hvl_ccb.dev(cube).switches

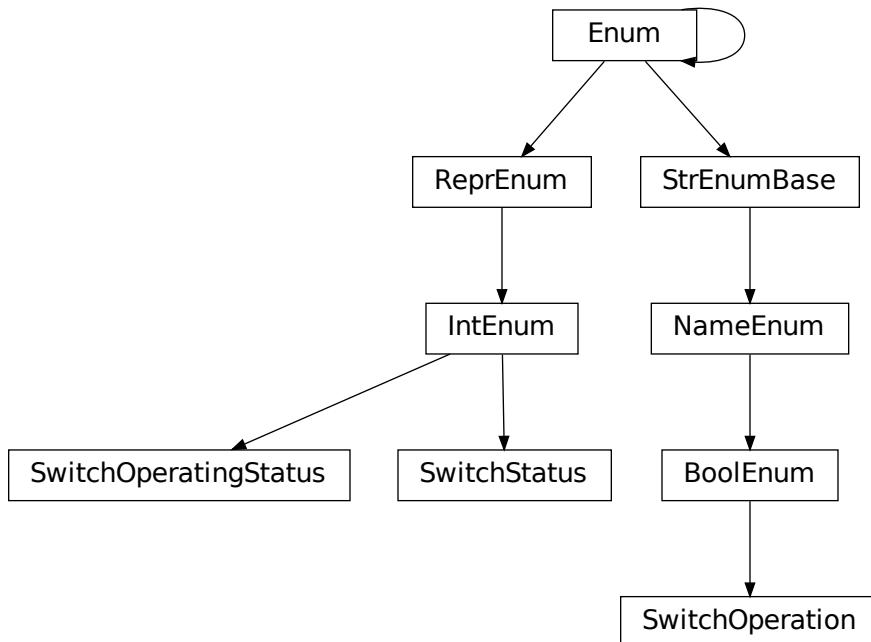

Switches of the different “Cubes”.

class SwitchOperatingStatus(*value=<no_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: **IntEnum**

Operating Status of a switch. Switch can be in auto or manual mode.

```
AUTO = 0
```

```
MANUAL = 1
```

```
class SwitchOperation(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *BoolEnum*

Operation of a switch in manual operating mode. Can be closed or opened.

```
CLOSE = True
```

```
OPEN = False
```

```
class SwitchStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *IntEnum*

Status of a switch. These are the possible values in the status integer e.g. in `_Switch.status`.

```
CLOSED = 1
```

```
ERROR = 3
```

```
INACTIVE = 0
```

```
OPEN = 2
```


Module contents

Cube package with implementation for system versions from 2019 on (new concept with hard-PLC Siemens S7-1500 as CPU).

`hvl_ccb.dev.ea_psi9000`

Submodules

`hvl_ccb.dev.ea_psi9000.ea_psi9000`

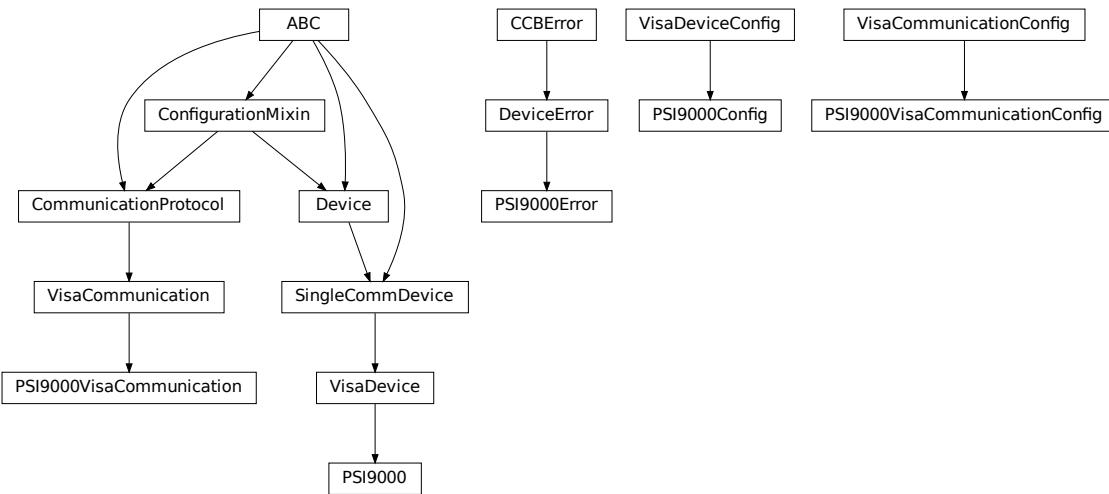

Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to now, all the testing was done with NI-Visa)

class `PSI9000`(*com*: `PSI9000VisaCommunication` | `PSI9000VisaCommunicationConfig` | *dict*, *dev_config*: `PSI9000Config` | *dict* | `None` = *None*)

Bases: `VisaDevice`

Elektro Automatik PSI 9000 power supply.

`MS_NOMINAL_CURRENT` = 2040

`MS_NOMINAL_VOLTAGE` = 80

`SHUTDOWN_CURRENT_LIMIT` = 0.1

`SHUTDOWN_VOLTAGE_LIMIT` = 0.1

`check_master_slave_config()` → *None*

Checks if the master / slave configuration and initializes if successful

Raises

`PSI9000Error` – if master-slave configuration failed

`static config_cls()`

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Return the default communication protocol for this device type, which is VisaCommunication.

Returns

the VisaCommunication class

get_output() → bool

Reads the current state of the DC output of the source. Returns True, if it is enabled, false otherwise.

Returns

the state of the DC output

get_system_lock() → bool

Get the current lock state of the system. The lock state is true, if the remote control is active and false, if not.

Returns

the current lock state of the device

get_ui_lower_limits() → tuple[float, float]

Get the lower voltage and current limits. A lower power limit does not exist.

Returns

Umin in V, Imin in A

get_ui_p_upper_limits() → tuple[float, float, float]

Get the upper voltage, current and power limits.

Returns

Umax in V, Imax in A, Pmax in W

get_voltage_current_setpoint() → tuple[float, float]

Get the voltage and current setpoint of the current source.

Returns

Uset in V, Iset in A

measure_voltage_current() → tuple[float, float]

Measure the DC output voltage and current

Returns

Umeas in V, Imeas in A

set_lower_limits(voltage_limit: float | None = None, current_limit: float | None = None) → None

Set the lower limits for voltage and current. After writing the values a check is performed if the values are set correctly.

Parameters

- **voltage_limit** – is the lower voltage limit in V
- **current_limit** – is the lower current limit in A

Raises

PSI9000Error – if the limits are out of range

set_output(target_onstate: bool) → None

Enables / disables the DC output.

Parameters

target_onstate – enable or disable the output power

Raises

PSI9000Error – if operation was not successful

set_system_lock(lock: bool) → None

Lock / unlock the device, after locking the control is limited to this class unlocking only possible when voltage and current are below the defined limits

Parameters

lock – True: locking, False: unlocking

set_upper_limits(voltage_limit: float | None = None, current_limit: float | None = None, power_limit: float | None = None) → None

Set the upper limits for voltage, current and power. After writing the values a check is performed if the values are set. If a parameter is left blank, the maximum configurable limit is set.

Parameters

- **voltage_limit** – is the voltage limit in V
- **current_limit** – is the current limit in A
- **power_limit** – is the power limit in W

Raises

PSI9000Error – if limits are out of range

set_voltage_current(volt: float, current: float) → None

Set voltage and current setpoints.

After setting voltage and current, a check is performed if writing was successful.

Parameters

- **volt** – is the setpoint voltage: 0.81.6 V (1.02 * 0-80 V) (absolute max, can be smaller if limits are set)
- **current** – is the setpoint current: 0..2080.8 A (1.02 * 0 - 2040 A) (absolute max, can be smaller if limits are set)

Raises

PSI9000Error – if the desired setpoint is out of limits

start() → None

Start this device.

stop() → None

Stop this device. Turns off output and lock, if enabled.

```
class PSI9000Config(spoll_interval: int | float = 0.5, spoll_start_delay: int | float = 2, power_limit: int | float = 43500, voltage_lower_limit: int | float = 0.0, voltage_upper_limit: int | float = 10.0, current_lower_limit: int | float = 0.0, current_upper_limit: int | float = 2040.0, wait_sec_system_lock: int | float = 0.5, wait_sec_settings_effect: int | float = 1, wait_sec_initialisation: int | float = 2)
```

Bases: *VisaDeviceConfig*

Elektro Automatik PSI 9000 power supply device class. The device is communicating over a VISA TCP socket.

Using this power supply, DC voltage and current can be supplied to a load with up to 2040 A and 80 V (using all four available units in parallel). The maximum power is limited by the grid, being at 43.5 kW available through the CEE63 power socket.

clean_values() → *None*

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

current_lower_limit: `int | float = 0.0`

Lower current limit in A, depending on the experimental setup.

current_upper_limit: `int | float = 2040.0`

Upper current limit in A, depending on the experimental setup.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

power_limit: `int | float = 43500`

Power limit in W depending on the experimental setup. With 3x63A, this is 43.5kW. Do not change this value, if you do not know what you are doing. There is no lower power limit.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

voltage_lower_limit: `int | float = 0.0`

Lower voltage limit in V, depending on the experimental setup.

voltage_upper_limit: `int | float = 10.0`

Upper voltage limit in V, depending on the experimental setup.

wait_sec_initialisation: `int | float = 2`**wait_sec_settings_effect:** `int | float = 1`**wait_sec_system_lock:** `int | float = 0.5`

exception PSI9000Error

Bases: *DeviceError*

Base error class regarding problems with the PSI 9000 supply.

class PSI9000VisaCommunication(*configuration*)

Bases: *VisaCommunication*

Communication protocol used with the PSI 9000 power supply.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

class PSI9000VisaCommunicationConfig(*host: str | IPv4Address | IPv6Address, interface_type: str | InterfaceType = InterfaceType.TCPIP_SOCKET, board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = ''*)

Bases: *VisaCommunicationConfig*

Visa communication protocol config dataclass with specification for the PSI 9000 power supply.

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

interface_type: str | InterfaceType = 1

Interface type of the VISA connection, being one of *InterfaceType*.

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

Module contents


Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to now, all the testing was done with NI-Visa)

`hvl_ccb.dev.fluke884x`

Submodules

`hvl_ccb.dev.fluke884x.base`

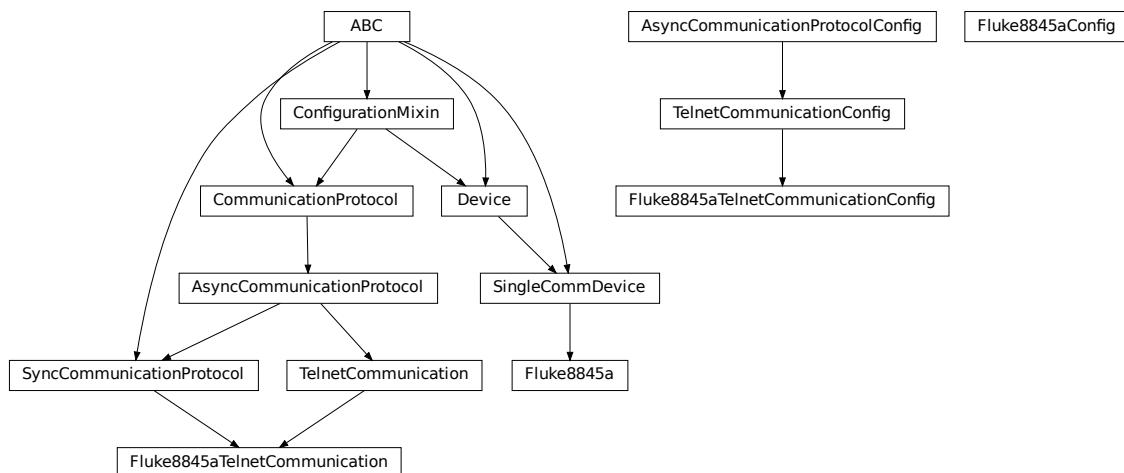

Python module for the Fluke8845a Multimeter. The communication to the device is through Telnet. 8845A/8846A Programmers Manual is available in the following link. All page numbers mentioned in this script refer to this manual. https://download.flukecal.com/pub/literature/8845A_pmeng0300.pdf

```
class Fluke8845a(com, dev_config=None)
```

Bases: `SingleCommDevice`

Device class to control Fluke8845a

`DISPLAY_MAX_LENGTH = 12`

`ac_current_range`

`ac_voltage_range`

`activate_remote_mode() → None`

Page 66

Places the Meter in the remote mode for RS-232 or Ethernet remote control. All front-panel keys, except the local key, are disabled.

clear_display_message() → *None*

Page 59

Clears the displayed message on the Meter's display.

clear_error_queue() → *None*

Page 62

Sets all bits to zero in the Meter's status byte register and all event registers. Also clears the error queue

static config_cls() → type[*Fluke8845aConfig*]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

current_filter

dc_current_range

dc_voltage_range

static default_com_cls() → type[*Fluke8845aTelnetCommunication*]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

property display_enable: bool

Page 59

get if the display is enabled or not fluke answer string “1” for ON and “0” for off bool(int(“1”)) = 1 and
bool(int(“0”)) = 0

Returns

bool enabled = True, else False

property display_message: str

Page 59

Retrieves the text sent to the Meter's display.

fetch() → float

Page 36

Transfer stored readings to output buffer

four_wire_resistance_range

frequency_aperture

property identification: str

Page 60

Queries “*IDN?” and returns the identification string of the connected device.

Returns

the identification string of the connected device e.g. “FLUKE, 8845A, 2540017, 08/02/10-11:53”

initiate_trigger() → *None*

Set trigger system to wait-for-trigger

measure() → float

Page 42

Taking measurement

Once the Meter has been configured for a measurement, the INITiate command causes the Meter to take a measurement when the trigger condition have been met. To process readings from the Meter's internal memory to the output buffer, send the Meter a FETCh? command.

property measurement_function: MeasurementFunction

input_function getter, query what the input function is

Raises

Fluke8845aUnknownCommandError – if the input function is unknown

period_aperture

reset() → *None*

Page 60

resets the meter to its power-up configuration

start() → *None*

Start this device as recommended by the manual

stop() → *None*

Stop this device. Disables access and closes the communication protocol.

trigger() → *None*

Causes the meter to trigger a measurement when paused

property trigger_delay: int

input_trigger_delay getter, query what the input trigger delay is in second answer format from Fluke: string, '+1.0000000E+00', so convert to float and then to int

Returns

input trigger delay in second

property trigger_source: TriggerSource

input_trigger_source getter, query what the input trigger source is

Raises

Fluke8845aUnknownCommandError – if the input trigger source is unknown

two_wire_resistance_range

voltage_filter

class Fluke8845aConfig(name: str = 'Fluke 1')

Bases: object

Config for Fluke8845a

name: the name of the device

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

name: str = 'Fluke 1'

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

class Fluke8845aTelnetCommunication(configuration)

Bases: *TelnetCommunication*, *SyncCommunicationProtocol*

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

query(*command: str*, *n_attempts_max: int | None = None*, *attempt_interval_sec: int | float | None = None*)

→ str

Send a command to the interface and handle the status message. Eventually raises an error.

Parameters

- **command** – Command to send
- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read as answer
- **attempt_interval_sec** – time between the reading attempts

Raises

Fluke8845aError – if the connection is broken

Returns

Answer from the interface

```
class Fluke8845aTelnetCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: int | float = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, host: str | ipaddress.IPv4Address | ipaddress.IPv6Address | NoneType = None, port: int = 3490, timeout: int | float = 0.2)
```

Bases: *TelnetCommunicationConfig*

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

port: int = 3490

Port at which Fluke 8845a is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns

a list of strings containing all required keys.

terminator: bytes = b'\r'

The terminator is CR

`hvl_ccb.dev.fluke884x.constants`

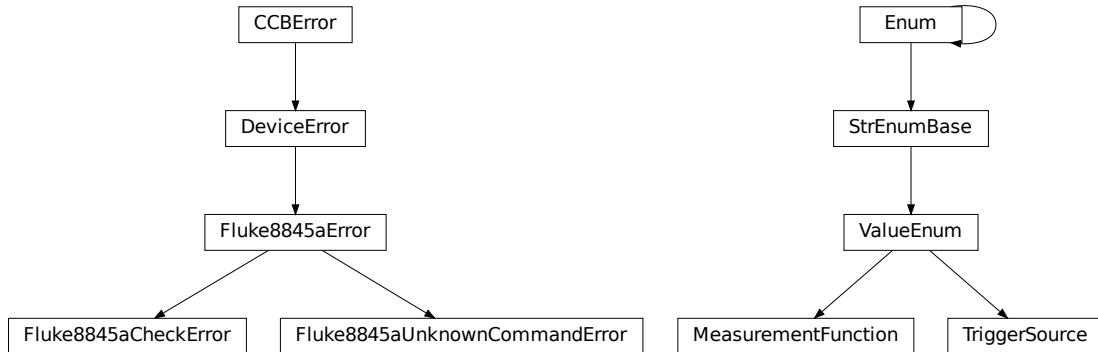

Constants, ValueEnum: MeasurementFunction and TriggerSoruce Descriptors for range, filter and aperture

exception Fluke8845aCheckError

Bases: *Fluke8845aError*

exception Fluke8845aError

Bases: *DeviceError*

exception Fluke8845aUnknownCommandError

Bases: *Fluke8845aError*

class MeasurementFunction(*value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None*)

Bases: *ValueEnum*

Page 40

Sets the Meter function. This command must be followed by the INIT and FETCh? commands to cause the meter to take a measurement.

CURRENT_AC = 'CURR:AC'

CURRENT_DC = 'CURR'

DIODE = 'DIOD'

FOUR_WIRE_RESISTANCE = 'FRES'

FREQUENCY = 'FREQ'

PERIOD = 'PER'

TWO_WIRE_RESISTANCE = 'RES'

VOLTAGE_AC = 'VOLT:AC'

VOLTAGE_DC = 'VOLT'

```
class TriggerSource(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *ValueEnum*

Page 57


BUS: Sets the Meter to expect a trigger through the IEEE-488 bus or upon execution of a *TRG command
IMM: Selects Meter's internal triggering system
EXT: Sets the Meter to sense triggers through the trigger jack on the rear panel of the Meter

```
BUS = 'BUS'
```

```
EXTERNAL = 'EXT'
```

```
IMMEDIATE = 'IMM'
```

hvl_ccb.dev.fluke884x.ranges

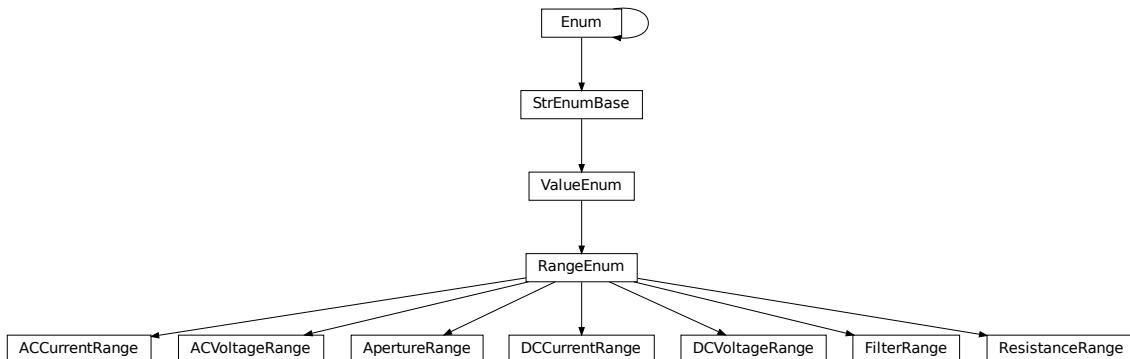

Ranges, RangeEnum for Fluke8845a device

```
class ACCurrentRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for AC current with unit Ampere

```
FOUR_HUNDRED_MILLI_AMPERE = 0.4
```

```
ONE_AMPERE = 1.0
```

```
ONE_HUNDRED_MILLI_AMPERE = 0.1
```

```
TEN_AMPERE = 10.0
```

```
TEN_MILLI_AMPERE = 0.01
```

```
THREE_AMPERE = 3.0
```

```
class ACVoltageRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for AC voltage with unit volt

HUNDRED_VOLT = 100.0

ONE_HUNDRED_MILLI_VOLT = 0.1

ONE_VOLT = 1.0

SEVEN_HUNDRED_FIFTY_VOLT = 750.0

TEN_VOLT = 10.0

```
class ApertureRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *RangeEnum*

Page 46

Sets the gate time for the frequency/period function to the value
10ms = 4 1/2 digits
100ms = 5 1/2 digits
1s = 6 1/2 digits

ONE_HUNDRED_MILLI_SECOND = 0.1

ONE_SECOND = 1.0

TEN_MILLI_SECOND = 0.01

```
class DCCurrentRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for DC current with unit Ampere

FOUR_HUNDRED_MILLI_AMPERE = 0.4

ONE_AMPERE = 1.0

ONE_HUNDRED_MICRO_AMPERE = 0.0001

ONE_HUNDRED_MILLI_AMPERE = 0.1

ONE_MILLI_AMPERE = 0.001

TEN_AMPERE = 10.0

TEN_MILLI_AMPERE = 0.01

THREE_AMPERE = 3.0

```
class DCVoltageRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for DC voltage with unit volt

```
HUNDRED_VOLT = 100.0
ONE_HUNDRED_MILLI_VOLT = 0.1
ONE_THOUSAND_VOLT = 1000.0
ONE_VOLT = 1.0
TEN_VOLT = 10.0

class FilterRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
```

Bases: *RangeEnum*

Page 47

```
Sets the appropriate filter for the frequency specified by <n>
High pass filter
For `VOLTAGE_AC`: <n> Hz to 300 kHz
For `CURRENT_AC`: <n> Hz to 10 kHz
parameters <n> = 3 slow filter
                  20 medium filter
                  200 fast filter
For `CURRENT_AC` and `VOLTAGE_AC`
```

```
FAST_FILTER = 200.0
```

```
MEDIUM_FILTER = 20.0
```

```
SLOW_FILTER = 3.0
```

```
class ResistanceRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                      boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for resistance with unit Ohm

```
ONE_HUNDRED_MILLION_OHM = 100000000.0
```

```
ONE_HUNDRED_OHM = 100.0
```

```
ONE_HUNDRED_THOUSAND_OHM = 100000.0
```

```
ONE_MILLION_OHM = 1000000.0
```

```
ONE_THOUSAND_OHM = 1000.0
```

```
TEN_MILLION_OHM = 10000000.0
```

```
TEN_THOUSAND_OHM = 10000.0
```


Module contents

Fluke 8845A multimeter implementation using Telnet communication

hvl_ccb.dev.fug

Submodules

hvl ccb.dev.fug.fug

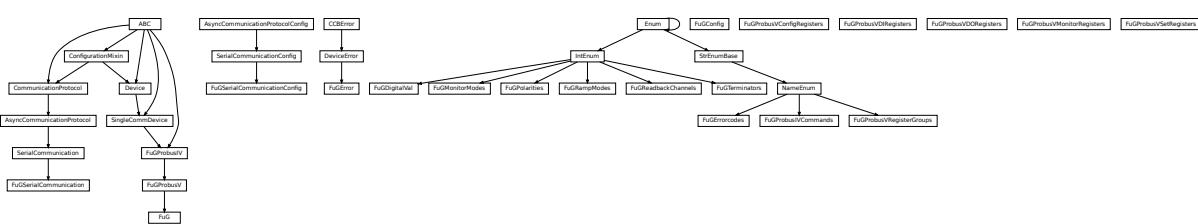

Device classes for “Probus V - ADDAT30” Interfaces which are used to control power supplies from FuG Elektronik GmbH

This interface is used for many FuG power units. Manufacturer homepage: <https://www.fug-elektronik.de>

The Professional Series of Power Supplies from FuG is a series of low, medium and high voltage direct current power supplies as well as capacitor chargers. The class FuG is tested with a HCK 800-20 000 in Standard Mode. The addressable mode is not implemented. Check the code carefully before using it with other devices. Manufacturer homepage: <https://www.fug-elektronik.de/netzgeraete/professional-series/>

The documentation of the interface from the manufacturer can be found here: https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus_V.zip

The provided classes support the basic and some advanced commands. The commands for calibrating the power supplies are not implemented, as they are only for very special porpoises and should not be used by "normal" customers.

```
class FuG(com, dev_config=None)
```

Bases: *FuGProbusV*

FuG power supply device class.

The power supply is controlled over a FuG ADDA Interface with the PROBUS V protocol

property config_status: *FuGProbusVConfigRegisters*

Returns the registers for the registers with the configuration and status values

Returns

FuGProbusVConfigRegisters

property current: *EuGProbusVSetRegisters*

Returns the registers for the current output.

Returns

property current monitor: *EuCProbusVMonitorRegisters*

Returns the registers for the current monitor

A typical usage will be “self.current_monitor.value” to measure the output current

Returns**property di: *FuGProbusVDIRegisters***

Returns the registers for the digital inputs

Returns

FuGProbusVDIRegisters

identify_device() → *None*

Identify the device nominal voltage and current based on its model number.

Raises

SerialCommunicationIOError – when communication port is not opened

property max_current: int | float

Returns the maximal current which could provided within the test setup

Returns**property max_current_hardware: int | float**

Returns the maximal current which could provided with the power supply

Returns**property max_voltage: int | float**

Returns the maximal voltage which could provided within the test setup

Returns**property max_voltage_hardware: int | float**

Returns the maximal voltage which could provided with the power supply

Returns**property on: *FuGProbusVDORegisters***

Returns the registers for the output switch to turn the output on or off

Returns

FuGProbusVDORegisters

property outX0: *FuGProbusVDORegisters*

Returns the registers for the digital output X0

Returns

FuGProbusVDORegisters

property outX1: *FuGProbusVDORegisters*

Returns the registers for the digital output X1

Returns

FuGProbusVDORegisters

property outX2: *FuGProbusVDORegisters*

Returns the registers for the digital output X2

Returns

FuGProbusVDORegisters

property outXCMD: *FuGProbusVDORegisters*

Returns the registers for the digital outputX-CMD

Returns

FuGProbusVDORegisters

start(*max_voltage=0, max_current=0*) → *None*

Opens the communication protocol and configures the device.

Parameters

- **max_voltage** – Configure here the maximal permissible voltage which is allowed in the given experimental setup
- **max_current** – Configure here the maximal permissible current which is allowed in the given experimental setup

property voltage: *FuGProbusVSetRegisters*

Returns the registers for the voltage output

Returns

property voltage_monitor: *FuGProbusVMonitorRegisters*

Returns the registers for the voltage monitor.

A typically usage will be “self.voltage_monitor.value” to measure the output voltage

Returns

class FuGConfig(*wait_sec_stop_commands: int | float = 0.5*)

Bases: *object*

Device configuration dataclass for FuG power supplies.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```

wait_sec_stop_commands: int | float = 0.5
    Time to wait after subsequent commands during stop (in seconds)

class FuGDigitalVal(value)
    Bases: IntEnum
    An enumeration.

    NO = 0
    OFF = 0
    ON = 1
    YES = 1

exception FuGError(*args, **kwargs)
    Bases: DeviceError
    Error with the FuG voltage source.

    errorcode: str
        Errorcode from the Probus, see documentation of Probus V chapter 5. Errors with three-digit errorcodes
        are thrown by this python module.

class FuGErrorcodes(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                     boundary=None)
    Bases: NameEnum

    The power supply can return an errorcode. These errorcodes are handled by this class. The original errorcodes
    from the source are with one or two digits, see documentation of Probus V chapter 5. All three-digit errorcodes
    are from this python module.

    E0 = ('no error', 'standard response on each command')
    E1 = ('no data available', 'Customer tried to read from GPIB but there were no data prepared. (IBIG50 sent command ~T2 to ADDA)')
    E10 = ('unknown SCPI command', 'This SCPI command is not implemented')
    E100 = ('Command is not implemented', 'You tried to execute a command, which is not implemented or does not exist')
    E106 = ('The rampstate is a read-only register', 'You tried to write data to the register, which can only give you the status of the ramping.')
    E11 = ('not allowed Trigger-on-Talk', 'Not allowed attempt to Trigger-on-Talk (~T1) while ADDA was in addressable mode.')
    E115 = ('The given index to select a digital value is out of range', 'Only integer values between 0 and 1 are allowed.')
    E12 = ('invalid argument in ~Tn command', 'Only ~T1 and ~T2 is implemented.')
    E125 = ('The given index to select a ramp mode is out of range', 'Only integer values between 0 and 4 are allowed.')
    E13 = ('invalid N-value', 'Register > K8 contained an invalid value. Error code is output on an attempt to query data with ? or ~T1')

```

```
E135 = ('The given index to select the readback channel is out of range', 'Only integer values between 0 and 6 are allowed.')

E14 = ('register is write only', 'Some registers can only be written to (i.e.> H0)')

E145 = ('The given value for the AD-conversion is unknown', 'Valid values for the ad-conversion are integer values from "0" to "7".')

E15 = ('string too long', 'i.e. serial number string too long during calibration')

E155 = ('The given value to select a polarity is out of range.', 'The value should be 0 or 1.')

E16 = ('wrong checksum', 'checksum over command string was not correct, refer also to 4.4 of the Probus V documentation')

E165 = ('The given index to select the terminator string is out of range', '')

E2 = ('unknown register type', "No valid register type after '>'")

E206 = ('This status register is read-only', 'You tried to write data to this register, which can only give you the actual status of the corresponding digital output.')

E306 = ('The monitor register is read-only', 'You tried to write data to a monitor, which can only give you measured data.')

E4 = ('invalid argument', 'The argument of the command was rejected .i.e. malformed number')

E5 = ('argument out of range', 'i.e. setvalue higher than type value')

E504 = ('Empty string as response', 'The connection is broken.')

E505 = ('The returned register is not the requested.', 'Maybe the connection is overburden.')

E6 = ('register is read only', 'Some registers can only be read but not written to. (i.e. monitor registers)')

E666 = ('You cannot overwrite the most recent error in the interface of the power supply. But, well: You created an error anyway...', '')

E7 = ('Receive Overflow', 'Command string was longer than 50 characters.')

E8 = ('EEPROM is write protected', 'Write attempt to calibration data while the write protection switch was set to write protected.')

E9 = ('address error', 'A non addressed command was sent to ADDA while it was in addressable mode (and vice versa).')

raise_()

class FuGMonitorModes(value)
Bases: IntEnum
An enumeration.
```

```

T1MS = 1
    15 bit + sign, 1 ms integration time

T200MS = 6
    typ. 19 bit + sign, 200 ms integration time

T20MS = 3
    17 bit + sign, 20 ms integration time

T256US = 0
    14 bit + sign, 256 us integration time

T40MS = 4
    17 bit + sign, 40 ms integration time

T4MS = 2
    15 bit + sign, 4 ms integration time

T800MS = 7
    typ. 20 bit + sign, 800 ms integration time

T80MS = 5
    typ. 18 bit + sign, 80 ms integration time

class FuGPolarities(value)
    Bases: IntEnum
    An enumeration.

    NEGATIVE = 1
    POSITIVE = 0

class FuGProbusIV(com, dev_config=None)
    Bases: SingleCommDevice, ABC
    FuG Probus IV device class

    Sends basic SCPI commands and reads the answer. Only the special commands and PROBUS IV instruction set is implemented.

    command(command: FuGProbusIVCommands, value=None) → str

        Parameters
            • command – one of the commands given within FuGProbusIVCommands
            • value – an optional value, depending on the command

        Returns
            a String if a query was performed

    static config_cls()
        Return the default configdataclass class.

        Returns
            a reference to the default configdataclass class

```

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

output_off() → None

Switch DC voltage output off.

reset() → None

Reset of the interface: All setvalues are set to zero

abstract start()

Open the associated communication protocol.

stop() → None

Close the associated communication protocol.

class FuGProbusIVCommands(*value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None*)

Bases: *NameEnum*

An enumeration.

ADMODE = ('S', (<enum 'FuGMonitorModes'>, <class 'int'>))

CURRENT = ('I', (<class 'int'>, <class 'float'>))

EXECUTE = ('X', None)

EXECUTEONX = ('G', (<enum 'FuGDigitalVal'>, <class 'int'>))

Wait for “X” to execute pending commands

ID = ('*IDN?', None)

OUTPUT = ('F', (<enum 'FuGDigitalVal'>, <class 'int'>))

POLARITY = ('P', (<enum 'FuGPolarities'>, <class 'int'>))

QUERY = ('?', None)

READBACKCHANNEL = ('N', (<enum 'FuGReadbackChannels'>, <class 'int'>))

RESET = ('=', None)

TERMINATOR = ('Y', (<enum 'FuGTerminators'>, <class 'int'>))

VOLTAGE = ('U', (<class 'int'>, <class 'float'>))

XOUTPUTS = ('R', <class 'int'>)

TODO: the possible values are limited to 0..13

class FuGProbusV(*com, dev_config=None*)

Bases: *FuGProbusIV*

FuG Probus V class which uses register based commands to control the power supplies

get_register(register: str) → str

get the value from a register

Parameters

register – the register from which the value is requested

Returns

the value of the register as a String

set_register(register: str, value: int | float | str) → None

generic method to set value to register

Parameters

- register – the name of the register to set the value
- value – which should be written to the register

class FuGProbusVConfigRegisters(fug, super_register: FuGProbusVRegisterGroups)

Bases: object

Configuration and Status values, acc. 4.2.5

property execute_on_x: FuGDigitalVal

status of Execute-on-X

Returns

FuGDigitalVal of the status

property most_recent_error: FuGErrorcodes

Reads the Error-Code of the most recent command

Return FuGError

Raises

FuGError – if code is not “E0”

property readback_data: FuGReadbackChannels

Preselection of readout data for Trigger-on-Talk

Returns

index for the readback channel

property srq_mask: int

SRQ-Mask, Service-Request Enable status bits for SRQ 0: no SRQ Bit 2: SRQ on change of status to CC
Bit 1: SRQ on change to CV

Returns

representative integer value

property srq_status: str

SRQ-Statusbyte output as a decimal number: Bit 2: PS is in CC mode Bit 1: PS is in CV mode

Returns

representative string

property status: str

Statusbyte as a string of 0/1. Combined status (compatibel to Probus IV), MSB first: Bit 7: I-REG Bit 6: V-REG Bit 5: ON-Status Bit 4: 3-Reg Bit 3: X-Stat (polarity) Bit 2: Cal-Mode Bit 1: unused Bit 0: SEL-D

Returns

string of 0/1

property terminator: *FuGTerminators*

Terminator character for answer strings from ADDA

Returns

FuGTerminators

class FuGProbusVDIRegisters(*fug, super_register:* FuGProbusVRegisterGroups)

Bases: object

Digital Inputs acc. 4.2.4

property analog_control: *FuGDigitalVal*

Returns

shows 1 if power supply is controlled by the analog interface

property calibration_mode: *FuGDigitalVal*

Returns

shows 1 if power supply is in calibration mode

property cc_mode: *FuGDigitalVal*

Returns

shows 1 if power supply is in CC mode

property cv_mode: *FuGDigitalVal*

Returns

shows 1 if power supply is in CV mode

property digital_control: *FuGDigitalVal*

Returns

shows 1 if power supply is digitally controlled

property on: *FuGDigitalVal*

Returns

shows 1 if power supply ON

property reg_3: *FuGDigitalVal*

For special applications.

Returns

input from bit 3-REG

property x_stat: *FuGPolarities*

Returns

polarity of HVPS with polarity reversal

class FuGProbusVDORegisters(*fug, super_register:* FuGProbusVRegisterGroups)

Bases: object

Digital outputs acc. 4.2.2

property out: int | *FuGDigitalVal*

Status of the output according to the last setting. This can differ from the actual state if output should only pulse.

Returns

FuGDigitalVal

property status: *FuGDigitalVal*

Returns the actual value of output. This can differ from the set value if pulse function is used.

Returns

FuGDigitalVal

class *FuGProbusVMonitorRegisters*(*fug, super_register: FuGProbusVRegisterGroups*)

Bases: *object*

Analog monitors acc. 4.2.3

property *adc_mode: FuGMonitorModes*

The programmed resolution and integration time of the AD converter

Returns

FuGMonitorModes

property *value: float*

Value from the monitor.

Returns

a float value in V or A

property *value_raw: float*

uncalibrated raw value from AD converter

Returns

float value from ADC

class *FuGProbusVRegisterGroups*(*value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None*)

Bases: *NameEnum*

An enumeration.

CONFIG = 'K'

INPUT = 'D'

MONITOR_I = 'M1'

MONITOR_V = 'M0'

OUTPUTONCMD = 'BON'

OUTPUTX0 = 'B0'

OUTPUTX1 = 'B1'

OUTPUTX2 = 'B2'

OUTPUTXCMD = 'BX'

SETCURRENT = 'S1'

SETVOLTAGE = 'S0'

class *FuGProbusVSetRegisters*(*fug, super_register: FuGProbusVRegisterGroups*)

Bases: *object*

Setvalue control acc. 4.2.1 for the voltage and the current output

property actualsetvalue: float

The actual valid set value, which depends on the ramp function.

Returns

actual valid set value

property high_resolution: FuGDigitalVal

Status of the high resolution mode of the output.

Return 0

normal operation

Return 1

High Res. Mode

property rampmode: FuGRampModes

The set ramp mode to control the setvalue.

Returns

the mode of the ramp as instance of FuGRampModes

property ramprate: float

The set ramp rate in V/s.

Returns

ramp rate in V/s

property rampstate: FuGDigitalVal

Status of ramp function.

Return 0

if final setvalue is reached

Return 1

if still ramping up

property setvalue: float

For the voltage or current output this setvalue was programmed.

Returns

the programmed setvalue

class FuGRampModes(value)

Bases: IntEnum

An enumeration.

FOLLOWRAMP = 1

Follow the ramp up- and downwards

IMMEDIATELY = 0

Standard mode: no ramp

ONLYUPWARDSOFFTOZERO = 4

Follow the ramp up- and downwards, if output is OFF set value is zero

RAMPUPWARDS = 2

Follow the ramp only upwards, downwards immediately

SPECIALRAMPUPWARDS = 3

Follow a special ramp function only upwards

```
class FuGReadbackChannels(value)
```

Bases: `IntEnum`

An enumeration.

CURRENT = 1

FIRMWARE = 5

RATEDCURRENT = 4

RATEDVOLTAGE = 3

SN = 6

STATUSBYTE = 2

VOLTAGE = 0

```
class FuGSerialCommunication(configuration)
```

Bases: `SerialCommunication`

Specific communication protocol implementation for FuG power supplies. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

query(*command*: str) → str

Send a command to the interface and handle the status message. Raises an error, if the answer starts with “E”.

Parameters

command – Command to send

Raises

`FuGError` – if the connection is broken or the error from the power source itself

Returns

Answer from the interface or empty string

```
class FuGSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',  
    encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:  
    Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int =  
    10, port: Optional[str] = None, baudrate: int = 9600, parity:  
    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =  
    <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,  
    hvl_ccb.comm.serial.SerialCommunicationStopbits] =  
    <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,  
    hvl_ccb.comm.serial.SerialCommunicationBytesize] =  
    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int,  
    float] = 3)
```

Bases: `SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for FuG power supplies is 9600 baud

bytesize: int | *SerialCommunicationBytesize* = 8

One byte is eight bits long

default_n_attempts_read_text_nonempty: int = 10

default number of attempts to read a non-empty text

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | *SerialCommunicationParity* = 'N'

FuG does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

stopbits: int | *SerialCommunicationStopbits* = 1

FuG uses one stop bit

terminator: bytes = b'\n'

The terminator is LF

timeout: int | float = 3

use 3 seconds timeout as default

wait_sec_read_text_nonempty: int | float = 0.5

default time to wait between attempts of reading a non-empty text

class FuGTerminators(*value*)

Bases: IntEnum

An enumeration.

CR = 3

```
CRLF = 0  
LF = 2  
LFCR = 1
```

Module contents

Device classes for “Probus V - ADDAT30” Interfaces which are used to control power supplies from FuG Elektronik GmbH

This interface is used for many FuG power units. Manufacturer homepage: <https://www.fug-elektronik.de>

The Professional Series of Power Supplies from FuG is a series of low, medium and high voltage direct current power supplies as well as capacitor chargers. The class FuG is tested with a HCK 800-20 000 in Standard Mode. The addressable mode is not implemented. Check the code carefully before using it with other devices. Manufacturer homepage: <https://www.fug-elektronik.de/netzgeraete/professional-series/>


The documentation of the interface from the manufacturer can be found here: https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus_V.zip

The provided classes support the basic and some advanced commands. The commands for calibrating the power supplies are not implemented, as they are only for very special porpoises and should not be used by “normal” customers.

hvl_ccb.dev.heinzinger

Submodules

`hvl_ccb.dev.heinzinger.base`

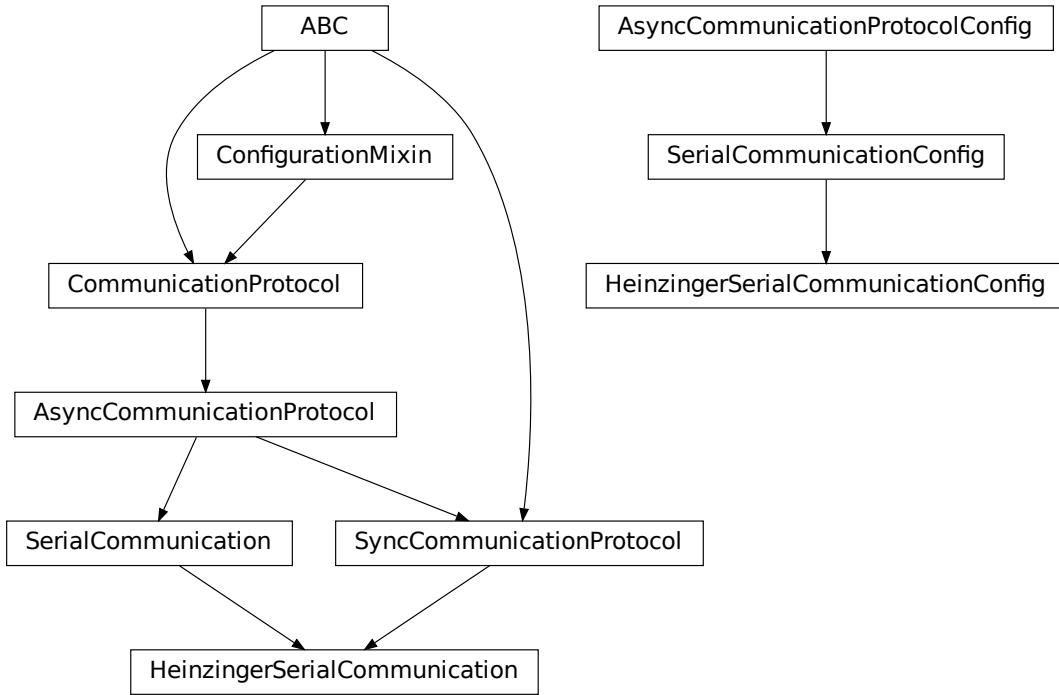

Base classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Interface Manual: https://www.heinzinger.com/assets/uploads/downloads/Handbuch_DigitalInterface_2021-12-14-V1.6.pdf

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class HeinzingerPNC is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCcap. Manufacturer homepage: <https://www.heinzinger.com/en/products/pnc-serie>

class HeinzingerSerialCommunication(configuration)

Bases: *SerialCommunication*, *SyncCommunicationProtocol*

Specific communication protocol implementation for Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class HeinzingerSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
    encoding_error_handling: str = 'strict',
    wait_sec_read_text_nonempty: Union[int, float] = 0.5,
    default_n_attempts_read_text_nonempty: int = 40, port:
    Optional[str] = None, baudrate: int = 9600, parity:
    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
    <SerialCommunicationParity.NONE: 'N'>, stopbits:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationStopbits] =
    <SerialCommunicationStopbits.ONE: 1>, bytesize:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationBytesize] =
    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
    Union[int, float] = 3)
```

Bases: *SerialCommunicationConfig*

baudrate: int = 9600

Baudrate for Heinzinger power supplies is 9600 baud

bytesize: int | SerialCommunicationBytesize = 8

One byte is eight bits long

default_n_attempts_read_text_nonempty: int = 40

increased to 40 default number of attempts to read a non-empty text

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

Heinzinger does not use parity

classmethod required_keys() → Sequence[str]


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
stopbits: int | SerialCommunicationStopbits = 1
    Heinzinger uses one stop bit
terminator: bytes = b'\n'
    The terminator is LF
timeout: int | float = 3
    use 3 seconds timeout as default
wait_sec_read_text_nonempty: int | float = 0.5
    default time to wait between attempts of reading a non-empty text
```

`hvl_ccb.dev.heinzinger.constants`

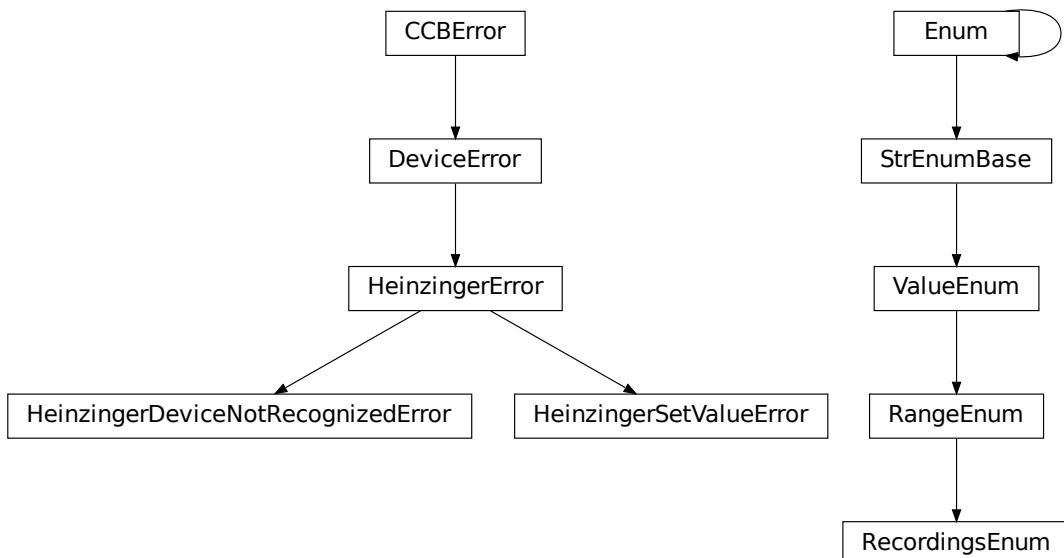

Constants, Heinzinger Digital Interface I/II and Heinzinger PNC power supply. Descriptors for errors

`exception HeinzingerDeviceNotRecognizedError`

Bases: `HeinzingerError`

Error indicating that the serial number of the device is not recognized.

`exception HeinzingerError`

Bases: `DeviceError`

General error with the Heinzinger PNC voltage source.

`exception HeinzingerSetValueError`

Bases: `HeinzingerError`

Error indicating that the value (current, voltage, ...) is not set correctly

```
class RecordingsEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *RangeEnum*

An enumeration.

EIGHT = 8.0


FOUR = 4.0

ONE = 1.0

SIXTEEN = 16.0

TWO = 2.0

hvl_ccb.dev.heinzinger.device

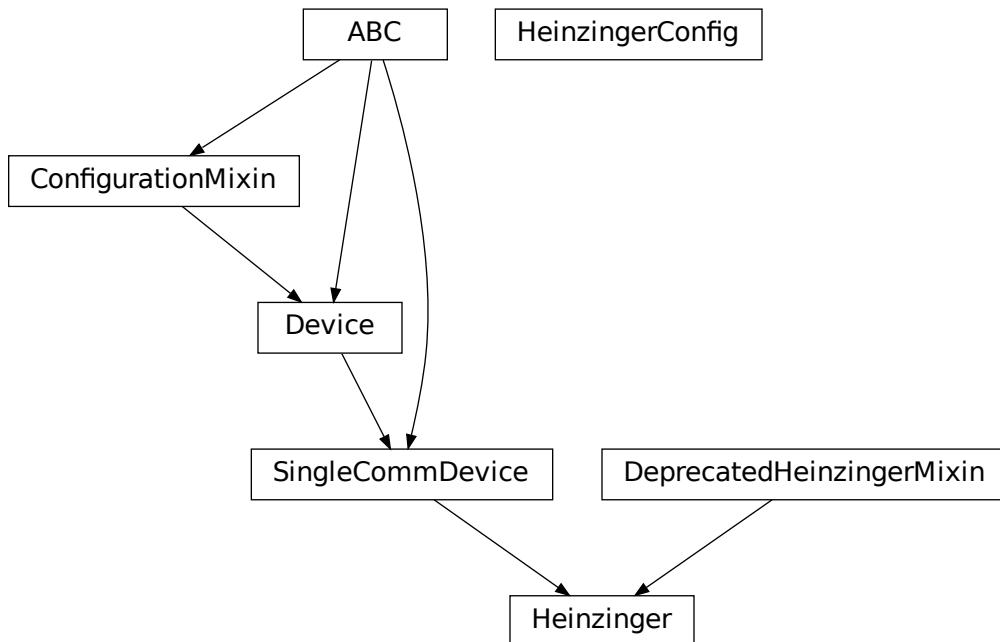

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Interface Manual: https://www.heinzinger.com/assets/uploads/downloads/Handbuch_DigitalInterface_2021-12-14-V1.6.pdf

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class Heinzinger is tested with different PNChp-types. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCcap. Manufacturer homepage: <https://www.heinzinger.com/en/products/pnc-serie>

class Heinzinger(*com, dev_config=None*)

Bases: *SingleCommDevice, DeprecatedHeinzingerMixin*

Heinzinger power supply device class. The power supply is controlled over a Heinzinger Digital Interface I/II

Sends basic SCPI commands and reads the answer. Only the standard instruction set from the manual is implemented.

static config_cls() → type[*HeinzingerConfig*]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property current: int | float

Ask the Device to measure its output current and return the measurement result in A.

Returns

measured current as float

Raises

SerialCommunicationIOError – when communication port is not opened

static default_com_cls() → type[*HeinzingerSerialCommunication*]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

get_interface_version() → str

Queries the version number of the digital interface.

Raises

SerialCommunicationIOError – when communication port is not opened

get_serial_number() → str

Ask the device for its serial number and returns the answer as a string.

Returns

string containing the device serial number

Raises

SerialCommunicationIOError – when communication port is not opened

identify_device() → *None*

Identify the device nominal voltage and current based on its serial number.

Raises

SerialCommunicationIOError – when communication port is not opened

property max_current: int | float

Maximal settable output current in A

property max_current.hardware: int | float

Maximal output current the hardware can provide in A

property max_voltage: int | float

Maximal settable output voltage in V

property max_voltage.hardware: int | float

Maximal output voltage the hardware can provide in V

property number_of_recordings: int

Queries the number of recordings the device is using for average value calculation.

Returns

int number of recordings

Raises

SerialCommunicationIOError – when communication port is not opened

property output: bool | None

Switch DC voltage output on and updates the output status.

Raises

SerialCommunicationIOError – when communication port is not opened

reset_interface() → None

Reset of the digital interface; only Digital Interface I: Power supply is switched to the Local-Mode (Manual operation)

Raises

SerialCommunicationIOError – when communication port is not opened

property set_current: int | float

Queries the set current of the Heinzinger PNC (not the measured current!) in A.

Raises

SerialCommunicationIOError – when communication port is not opened

property set_voltage: int | float

Queries the set voltage of the Heinzinger PNC (not the measured voltage!) in V.

Raises

SerialCommunicationIOError – when communication port is not opened

start() → None

Opens the communication protocol and configures the device.

Raises

SerialCommunicationIOError – when communication port cannot be opened.

stop() → None

Stop the device. Closes also the communication protocol.

property voltage: int | float

Ask the Device to measure its output voltage and return the measurement result in V.

Returns

measured voltage as float in V

Raises

SerialCommunicationIOError – when communication port is not opened

class HeinzingerConfig(default_number_of_recordings: int | RecordingsEnum = 1, number_of_decimals: int = 6, wait_sec_stop_commands: int | float = 0.5)

Bases: object

Device configuration dataclass for Heinzinger power supplies.

clean_values()

default_number_of_recordings: int | RecordingsEnum = 1

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

number_of_decimals: int = 6**classmethod optional_defaults()** → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

wait_sec_stop_commands: int | float = 0.5

Time to wait after subsequent commands during stop (in seconds)

hvl_ccb.dev.heinzinger.mixin**DeprecatedHeinzingerMixin**

Mixin class for Heinzinger's deprecated methods.Raise

DeprecationWarning, will be removed in the next release

class DeprecatedHeinzingerMixin

Bases: object

static get_number_of_recordings() → *None***static measure_current()** → *None*

```
static measure_voltage() → None
static output_off() → None
static output_on() → None
static set_number_of_recordings() → None
```

Module contents

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.


The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Interface Manual: https://www.heinzinger.com/assets/uploads/downloads/Handbuch_DigitalInterface_2021-12-14-V1.6.pdf

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class HeinzingerPNC is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCcap. Manufacturer homepage: <https://www.heinzinger.com/en/products/pnc-serie>

hvl_ccb.dev.highland_t560

Submodules

hvl_ccb.dev.highland_t560.base

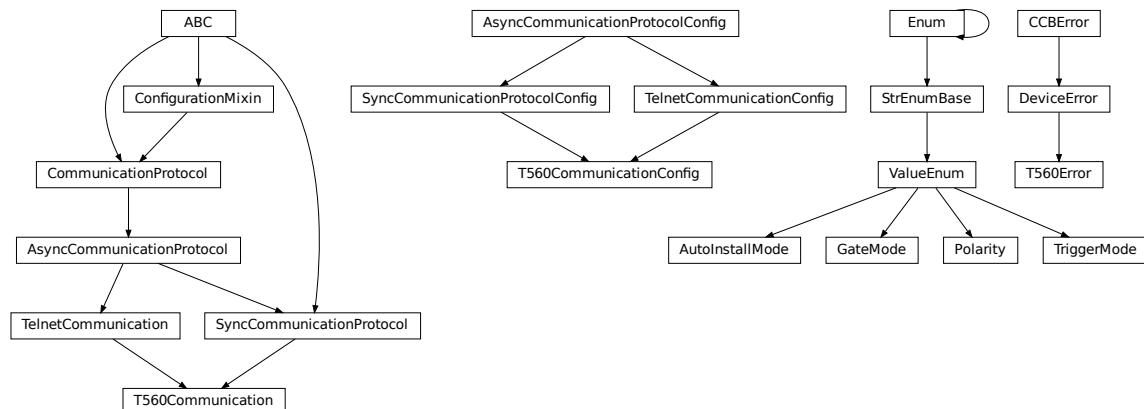

Module containing base device and communication classes and enums.

Communication with device is performed via its ethernet port and a Telnet connection.

```
class AutoInstallMode(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: `ValueEnum`

Modes for installing configuration settings to the device.

```
INSTALL = 1
```

```
OFF = 0
QUEUE = 2

class GateMode(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                boundary=None)
Bases: ValueEnum
Available T560 gate modes
INPUT = 'INP'
OFF = 'OFF'
OUTPUT = 'OUT'

class Polarity(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                boundary=None)
Bases: ValueEnum
Possible channel polarity states
ACTIVE_HIGH = 'POS'
ACTIVE_LOW = 'NEG'

class T560Communication(configuration)
Bases: SyncCommunicationProtocol, TelnetCommunication
Communication class for T560. It uses a TelnetCommunication with the SyncCommunicationProtocol
static config_cls()
Return the default configdataclass class.

    Returns
        a reference to the default configdataclass class

query(command: str, n_attempts_max: int | None = None, attempt_interval_sec: int | float | None = None)
    → str
Send a command to the device and handle the response.
For device setting queries, response will be ‘OK’ if successful, or ‘?’ if setting cannot be carried out,
raising an error.

    Parameters
        • command – Command string to be sent
        • n_attempts_max – Amount of attempts how often a non-empty text is tried to be read as
            answer
        • attempt_interval_sec – time between the reading attempts

    Raises
        T560Error – if no response is received, or if the device responds with an error message.

    Returns
        Response from the device.
```

```
class T560CommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: int | float = 0.5, default_n_attempts_read_text_nonempty: int = 10, host: str | ipaddress.IPv4Address | ipaddress.IPv6Address | NoneType = None, port: int = 2000, timeout: int | float = 0.2)
```

Bases: *SyncCommunicationProtocolConfig*, *TelnetCommunicationConfig*

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

port: int = 2000

Port at which the host is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

terminator: bytes = b'\r'

The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations. This defines the end of a single line.

exception T560Error

Bases: *DeviceError*

T560 related errors.

```
class TriggerMode(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *ValueEnum*

Available T560 trigger modes

COMMAND = 'REM'


EXT_FALLING_EDGE = 'NEG'

```
EXT_RISING_EDGE = 'POS'  
INT_SYNTHESIZER = 'SYN'  
OFF = 'OFF'
```

hvl_ccb.dev.highland_t560.channel

Module for controlling pulse output channels A, B, C and D.

hvl_ccb.dev.highland_t560.device

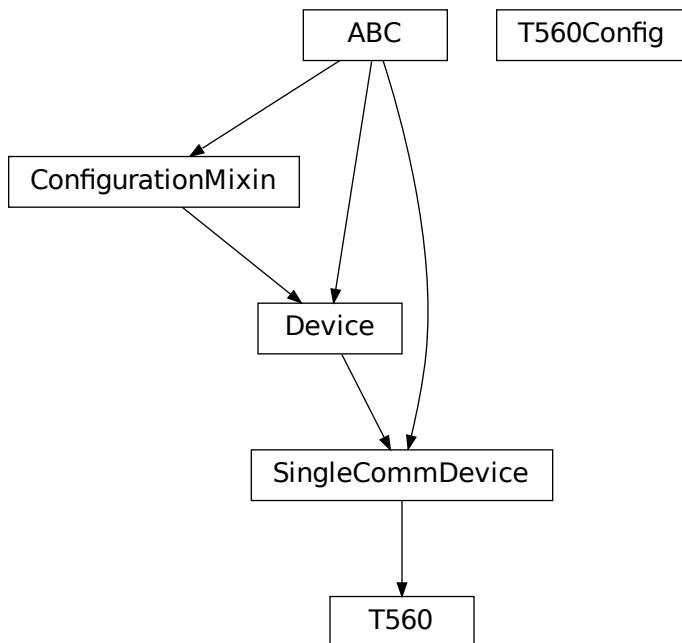

including TRIG, CLOCK and GATE I/Os.

Module for controlling device,

```
class T560(com, dev_config=None)
```

Bases: `SingleCommDevice`

activate_clock_output()

Outputs 10 MHz clock signal

property auto_install_mode: `AutoInstallMode`

Check the autoinstall settings of the T560. The autoinstall mode sets how changes to device settings are applied. See manual section 4.7.2 for more information about these modes.

property ch_a: `_Channel`

Channel A of T560

```
property ch_b: _Channel
    Channel B of T560

property ch_c: _Channel
    Channel C of T560

property ch_d: _Channel
    Channel D of T560

static config_cls()
    Return the default configdataclass class.

Returns
    a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

Returns
    the type of the standard communication protocol for this device

disarm_trigger()
    Disarm DDG by disabling all trigger sources.

fire_trigger()
    Fire a software trigger.

property frequency: float
    The frequency of the timing cycle in Hz.

property gate_mode: GateMode
    Check the mode setting of the GATE I/O port.

property gate_polarity: Polarity
    Check the polarity setting of the GATE I/O port.

load_device_configuration()
    Load the settings saved in nonvolatile memory.

property period: float
    The period of the timing cycle (time between triggers) in seconds.

save_device_configuration()
    Save the current settings to nonvolatile memory.

property trigger_level
    Get external trigger level.

property trigger_mode
    Get device trigger source.

use_external_clock()
    Finds and accepts an external clock signal to the CLOCK input

class T560Config
    Bases: object
        auto_install_mode = 1
```

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.


Module contents

This module establishes methods for interfacing with the Highland Technology T560-2 via its ethernet adapter with a telnet communication protocol.

The T560 is a small digital delay & pulse generator. It outputs up to four individually timed pulses with 10-ps precision, given an internal or external trigger.

This module introduces methods for configuring channels, gating, and triggering. Further documentation and a more extensive command list may be obtained from:

<https://www.highlandtechnology.com/DSS/T560DS.shtml>

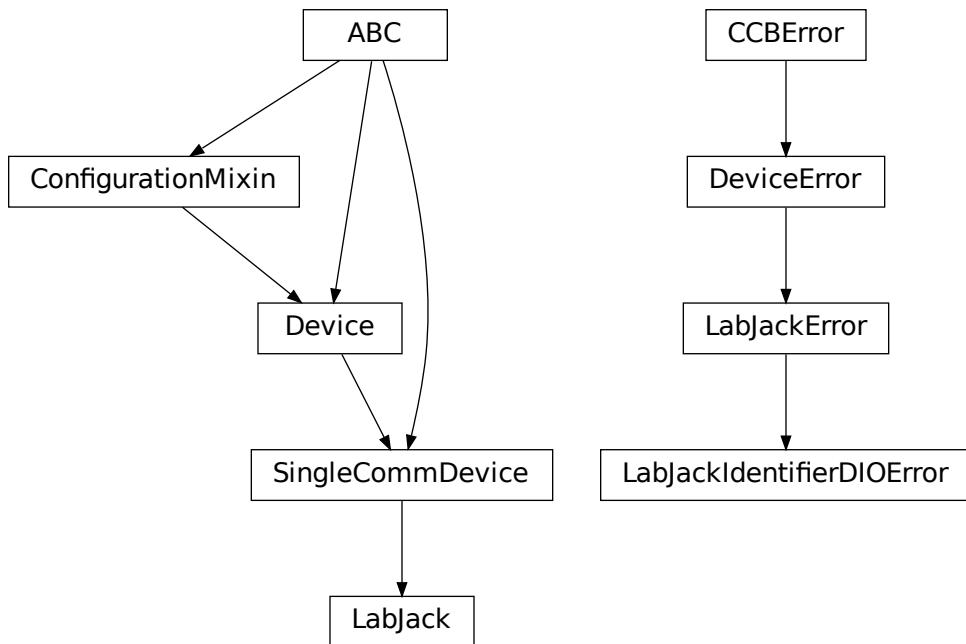hvl_ccb.dev.labjack**Submodules****hvl_ccb.dev.labjack.labjack**

A
LabJack T-series devices wrapper around the LabJack's LJM Library; see <https://labjack.com/ljm> . The wrapper was originally developed and tested for a LabJack T7-PRO device.

Extra installation

To use this LabJack T-series devices wrapper:

1. install the hvl_ccb package with a labjack extra feature:

```
$ pip install "hvl_ccb[labjack]"
```

this will install the Python bindings for the library.

2. install the library - follow instruction in <https://labjack.com/support/software/installers/ljm> .

class **LabJack**(*com*, *dev_config=None*)

Bases: *SingleCommDevice*

LabJack Device.

This class is tested with a LabJack T7-Pro and should also work with T4 and T7 devices communicating through the LJM Library. Other or older hardware versions and variants of LabJack devices are not supported.

```
class AInRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: [StrEnumBase](#)

An enumeration.

ONE = 1.0

ONE_HUNDREDTH = 0.01

ONE_TENTH = 0.1

TEN = 10.0

property value: float

```
class BitLimit(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: [IntEnum](#)

Maximum integer values for clock settings

THIRTY_TWO_BIT = 4294967295

```
class CalMicroAmpere(value=<no_arg>, names=None, module=None, qualname=None, type=None,  
start=1, boundary=None)
```

Bases: [Enum](#)

Pre-defined microampere (uA) values for calibration current source query.

TEN = '10uA'

TWO_HUNDRED = '200uA'

```
class CjcType(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: [NameEnum](#)

CJC slope and offset

internal = (1, 0)

lm34 = (55.56, 255.37)

```
class ClockFrequency(value=<no_arg>, names=None, module=None, qualname=None, type=None,  
start=1, boundary=None)
```

Bases: [IntEnum](#)

Available clock frequencies, in Hz

FIVE_MHZ = 5000000

FORTY_MHZ = 40000000

MAXIMUM = 80000000

MINIMUM = 312500

```

TEN_MHZ = 10000000
TWELVE_HUNDRED_FIFTY_KHZ = 1250000
TWENTY_FIVE_HUNDRED_KHZ = 2500000
TWENTY_MHZ = 20000000

DIOChannel
alias of TSeriesDIOChannel

class DIOStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                 boundary=None)
Bases: IntEnum
State of a digital I/O channel.

HIGH = 1
LOW = 0

class DeviceType(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
Bases: AutoNumberNameEnum
LabJack device types.

Can be also looked up by ambiguous Product ID (p_id) or by instance name: `python
LabJackDeviceType(4) is LabJackDeviceType('T4')` 

ANY = 1
T4 = 2
T7 = 3
T7_PRO = 4

classmethod get_by_p_id(p_id: int) → DeviceType | list[DeviceType]
Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

Parameters
    p_id – Product ID of a LabJack device
Returns
    Instance or list of instances of LabJackDeviceType
Raises
    ValueError – when Product ID is unknown

class TemperatureUnit(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                      start=1, boundary=None)
Bases: NameEnum
Temperature unit (to be returned)

C = 1
F = 2
K = 0

```

```
class ThermocoupleType(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                       start=1, boundary=None)
```

Bases: [NameEnum](#)

Thermocouple type; NONE means disable thermocouple mode.

C = 30

E = 20

J = 21

K = 22

NONE = 0

PT100 = 40

PT1000 = 42

PT500 = 41

R = 23

S = 25

T = 24

```
config_high_pulse(address: str | TSeriesDIOChannel, t_start: int | float, t_width: int | float, n_pulses: int
                  = 1) → None
```

Configures one FIO channel to send a timed HIGH pulse. Configure multiple channels to send pulses with relative timing accuracy. Times have a maximum resolution of 1e-7 seconds @ 10 MHz. :param address: FIO channel: [T7] FIO0;2;3;4;5. [T4] FIO6;7. :raises LabJackError if address is not supported. :param t_start: pulse start time in seconds. :raises ValueError: if t_start is negative or would exceed the clock period. :param t_width: duration of high pulse, in seconds. :raises ValueError: if t_width is negative or would exceed the clock period. :param n_pulses: number of pulses to be sent; single pulse default. :raises TypeError if n_pulses is not of type int. :raises ValueError if n_pulses is negative or exceeds the 32bit limit.

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

```
disable_pulses(*addresses: str | TSeriesDIOChannel | None) → None
```

Disable previously configured pulse channels. :param addresses: tuple of FIO addresses. All channels disabled if no argument is passed.

```
enable_clock(clock_enabled: bool) → None
```

Enable/disable LabJack clock to configure or send pulses. :param clock_enabled: True -> enable, False -> disable. :raises TypeError: if clock_enabled is not of type bool

```
get_ain(*channels: int) → float | Sequence[float]
```

Read currently measured value (voltage, resistance, ...) from one or more of analog inputs.

Parameters

channels – AIN number or numbers (0..254)

Returns

the read value (voltage, resistance, ...) as `float` or `tuple` of them in case multiple channels given

get_cal_current_source(*name: str | CalMicroAmpere*) → `float`

This function will return the calibration of the chosen current source, this ist not a measurement!

The value was stored during fabrication.

Parameters

name – ‘200uA’ or ‘10uA’ current source

Returns

calibration of the chosen current source in ampere

get_clock() → `dict[str, int | float]`

Return clock settings read from LabJack.

get_digital_input(*address: str | TSeriesDIOChannel*) → `DIOStatus`

Get the value of a digital input.

allowed names for T7 (Pro): FIO0 - FIO7, EIO0 - EIO 7, CIO0- CIO3, MIO0 - MIO2 :param address: name of the output -> ‘FIO0’ :return: HIGH when *address* DIO is high, and LOW when *address* DIO is low

get_product_id() → `int`

This function returns the product ID reported by the connected device.

Attention: returns 7 for both T7 and T7-Pro devices!

Returns

integer product ID of the device

get_product_name(*force_query_id=False*) → `str`

This function will return the product name based on product ID reported by the device.

Attention: returns “T7” for both T7 and T7-Pro devices!

Parameters

force_query_id – boolean flag to force `get_product_id` query to device instead of using cached device type from previous queries.

Returns

device name string, compatible with `LabJack.DeviceType`

get_product_type(*force_query_id: bool = False*) → `DeviceType`

This function will return the device type based on reported device type and in case of unambiguity based on configuration of device’s communication protocol (e.g. for “T7” and “T7_PRO” devices), or, if not available first matching.

Parameters

force_query_id – boolean flag to force `get_product_id` query to device instead of using cached device type from previous queries.

Returns

`DeviceType` instance

Raises

`LabJackIdentifierDIOError` – when read Product ID is unknown

get_sbus_rh(*number: int*) → float

Read the relative humidity value from a serial SBUS sensor.

Parameters

number – port number (0..22)

Returns

relative humidity in %RH

get_sbus_temp(*number: int*) → float

Read the temperature value from a serial SBUS sensor.

Parameters

number – port number (0..22)

Returns

temperature in Kelvin

get_serial_number() → int

Returns the serial number of the connected LabJack.

Returns

Serial number.

read_resistance(*channel: int*) → float

Read resistance from specified channel.

Parameters

channel – channel with resistor

Returns

resistance value with 2 decimal places

read_thermocouple(*pos_channel: int*) → float

Read the temperature of a connected thermocouple.

Parameters

pos_channel – is the AIN number of the positive pin

Returns

temperature in specified unit

send_pulses(**addresses: str | TSeriesDIOChannel*) → *None*

Sends pre-configured pulses for specified addresses. :param *addresses*: tuple of FIO addresses :raises *LabJackError* if an address has not been configured.

set_ain_differential(*pos_channel: int, differential: bool*) → *None*

Sets an analog input to differential mode or not. T7-specific: For base differential channels, positive must be even channel from 0-12 and negative must be positive+1. For extended channels 16-127, see Mux80 datasheet.

Parameters

- **pos_channel** – is the AIN number (0..12)
- **differential** – True or False

Raises

LabJackError – if parameters are unsupported

set_ain_range(channel: int, vrage: Real | AInRange) → None

Set the range of an analog input port.

Parameters

- **channel** – is the AIN number (0..254)
- **vrage** – is the voltage range to be set

set_ain_resistance(channel: int, vrage: Real | AInRange, resolution: int) → None

Set the specified channel to resistance mode. It utilized the 200uA current source of the LabJack.

Parameters

- **channel** – channel that should measure the resistance
- **vrage** – voltage range of the channel
- **resolution** – resolution index of the channel T4: 0-5, T7: 0-8, T7-Pro 0-12

set_ain_resolution(channel: int, resolution: int) → None

Set the resolution index of an analog input port.

Parameters

- **channel** – is the AIN number (0..254)
- **resolution** – is the resolution index within 0...`get_product_type().ain_max_resolution` range; 0 will set the resolution index to default value.

set_ain_thermocouple(pos_channel: int, thermocouple: None | str | ThermocoupleType, cjc_address: int = 60050, cjc_type: str | CjcType = CjcType.internal, vrage: Real | AInRange = AInRange.ONE_HUNDREDTH, resolution: int = 10, unit: str | TemperatureUnit = TemperatureUnit.K) → None

Set the analog input channel to thermocouple mode.

Parameters

- **pos_channel** – is the analog input channel of the positive part of the differential pair
- **thermocouple** – None to disable thermocouple mode, or string specifying the thermocouple type
- **cjc_address** – modbus register address to read the CJC temperature
- **cjc_type** – determines cjc slope and offset, ‘internal’ or ‘lm34’
- **vrage** – measurement voltage range
- **resolution** – resolution index (T7-Pro: 0-12)
- **unit** – is the temperature unit to be returned (‘K’, ‘C’ or ‘F’)

Raises

LabJackError – if parameters are unsupported

set_analog_output(channel: int, value: int | float) → None

Set the voltage of a analog output port

Parameters

- **channel** – DAC channel number 1/0
- **value** – The output voltage value 0-5 Volts int/float

set_clock(*clock_frequency*: Number | ClockFrequency = 10000000, *clock_period*: Number = 1) → *None*

Configure LabJack clock for pulse out feature. :param *clock_frequency*: clock frequency in Hz; default 10 MHz for base 10. :raises ValueError: if *clock_frequency* is not allowed (see ClockFrequency). :param *clock_period*: clock roll time in seconds; default 1s, 0 for max. :raises ValueError: if *clock_period* exceeds the 32bit tick limit. Clock period determines pulse spacing when using multi-pulse settings. Ensure period exceeds maximum intended pulse end time.

set_digital_output(*address*: str, *state*: int | DIOStatus) → *None*

Set the value of a digital output.

Parameters

- **address** – name of the output -> ‘FIO0’
- **state** – state of the output -> *DIOStatus* instance or corresponding *int* value

start() → *None*

Start the Device.

stop() → *None*

Stop the Device.

exception LabJackError

Bases: *DeviceError*

General Error for the LabJack device.

exception LabJackIdentifierDIOError

Bases: *LabJackError*

Error indicating a wrong DIO identifier

Module contents

A LabJack T-series devices wrapper around the LabJack’s LJM Library; see <https://labjack.com/ljm> . The wrapper was originally developed and tested for a LabJack T7-PRO device.

Extra installation

To use this LabJack T-series devices wrapper:

1. install the hvl_ccb package with a labjack extra feature:

```
$ pip install "hvl_ccb[labjack]"
```


this will install the Python bindings for the library.

2. install the library - follow instruction in <https://labjack.com/support/software/installers/ljm> .

hvl_ccb.dev.lauda

Submodules

hvl_ccb.dev.lauda.lauda

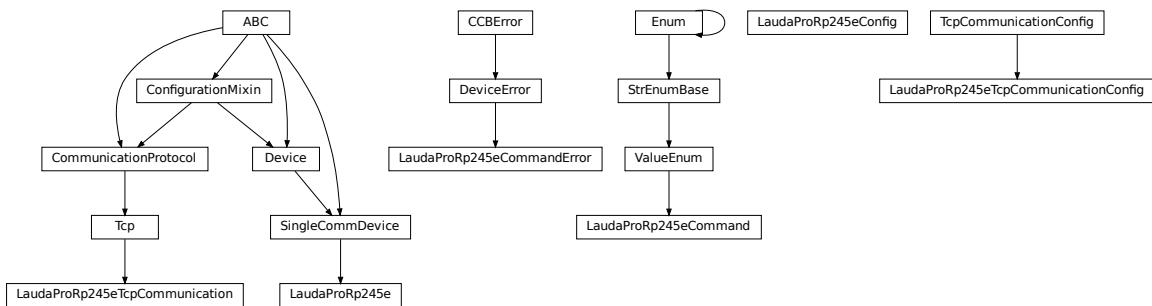

Device class for controlling a Lauda PRO RP245E, circulation chiller over TCP.

class LaudaProRp245e(com, dev_config=None)

Bases: *SingleCommDevice*

Lauda RP245E circulation chiller class.

static config_cls() → type[*LaudaProRp245eConfig*]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

continue_ramp() → str

Continue current ramp program.

Returns

reply of the device to the last call of “query”

static default_com_cls() → type[*LaudaProRp245eTcpCommunication*]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

get_bath_temp() → float

:return : float value of measured lauda bath temp in °C

get_device_type() → str

:return : Connected Lauda device type (for connection/com test)

pause() → str

Stop temperature control and pump.

Returns

reply of the device to the last call of “query”

pause_ramp() → str

Pause current ramp program.

Returns

reply of the device to the last call of “query”

reset_ramp() → str

Delete all segments from current ramp program.

Returns

reply of the device to the last call of “query”

run() → str

Start temperature control & pump.

Returns

reply of the device to the last call of “query”

set_control_mode(mod: int | ExtControlModeEnum = ExtControlModeEnum.INTERNAL) → str

Define control mode. 0 = INTERNAL (control bath temp), 1 = EXPT100 (pt100 attached to chiller), 2 = ANALOG, 3 = SERIAL, 4 = USB, 5 = ETH (to be used when passing the ext. temp. via ethernet) (temperature then needs to be passed every second, when not using options 3, 4, or 5)

Parameters

mod – temp control mode (control internal temp or external temp).

Returns

reply of the device to the last call of “query” (“OK”, if command was recognized”)

set_external_temp(external_temp: float = 20.0) → str

Pass value of external controlled temperature. Should be done every second, when control of external temperature is active. Has to be done right before control of external temperature is activated.

Parameters

external_temp – current value of external temperature to be controlled.

Returns

reply of the device to the last call of “query”

set_pump_level(pump_level: int = 6) → str

Set pump level Raises ValueError, if pump level is invalid.

Parameters

pump_level – pump level.

Returns

reply of the device to the last call of “query”

set_ramp_iterations(num: int = 1) → str

Define number of ramp program cycles.

Parameters

num – number of program cycles to be performed.

Returns

reply of the device to the last call of “query”

set_ramp_program(program: int = 1) → str

Define ramp program for following ramp commands. Raises ValueError if maximum number of ramp programs (5) is exceeded.

Parameters

program – Number of ramp program to be activated for following commands.

Returns

reply of the device to the last call of “query”

set_ramp_segment(*temp*: float = 20.0, *dur*: int = 0, *tol*: float = 0.0, *pump*: int = 6) → str

Define segment of current ramp program - will be attached to current program. Raises ValueError, if pump level is invalid.

Parameters

- **temp** – target temperature of current ramp segment
- **dur** – duration in minutes, in which target temperature should be reached
- **tol** – tolerance at which target temperature should be reached (for 0.00, next segment is started after dur has passed).
- **pump** – pump level to be used for this program segment.

Returns

reply of the device to the last call of “query”

set_temp_set_point(*temp_set_point*: float = 20.0) → str

Define temperature set point

Parameters

temp_set_point – temperature set point.

Returns

reply of the device to the last call of “query”

start() → *None*

Start this device.

start_ramp() → str

Start current ramp program.

Returns

reply of the device to the last call of “query”

stop() → *None*

Stop this device. Disables access and closes the communication protocol.

stop_ramp() → str

Stop current ramp program.

Returns

reply of the device to the last call of “query”

validate_pump_level(*level*: int)

Validates pump level. Raises ValueError, if pump level is incorrect. :param level: pump level, integer

class LaudaProRp245eCommand(*value*=<no_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *ValueEnum*

Commands for Lauda PRO RP245E Chiller Command strings most often need to be complimented with a parameter (attached as a string) before being sent to the device. Commands implemented as defined in “Lauda Betriebsanleitung fuer PRO Badthermostate und Umwaelzthermostate” pages 42 - 49

```
BATH_TEMP = 'IN_PV_00'  
    Request internal bath temperature  
COM_TIME_OUT = 'OUT_SP_08_'  
    Define communication time out  
CONT_MODE = 'OUT_MODE_01_'  
    Set control mode 1=internal, 2=ext. analog, 3=ext. serial, 4=USB, 5=ethernet  
DEVICE_TYPE = 'TYPE'  
    Request device type  
EXTERNAL_TEMP = 'OUT_PV_05_'  
    Pass on external controlled temperature  
LOWER_TEMP = 'OUT_SP_05_'  
    Define lower temp limit  
OPERATION_MODE = 'OUT_SP_02_'  
    Define operation mode  
PUMP_LEVEL = 'OUT_SP_01_'  
    Define pump level 1-8  
RAMP_CONTINUE = 'RMP_CONT'  
    Continue a paused ramp program  
RAMP_DELETE = 'RMP_RESET'  
    Reset a selected ramp program  
RAMP_ITERATIONS = 'RMP_OUT_02_'  
    Define how often a ramp program should be iterated  
RAMP_PAUSE = 'RMP_PAUSE'  
    Pause a selected ramp program  
RAMP_SELECT = 'RMP_SELECT_'  
    Select a ramp program (target for all further ramp commands)  
RAMP_SET = 'RMP_OUT_00_'  
    Define parameters of a selected ramp program  
RAMP_START = 'RMP_START'  
    Start a selected ramp program  
RAMP_STOP = 'RMP_STOP'  
    Stop a running ramp program  
START = 'START'  
    Start temp control (pump and heating/cooling)  
STOP = 'STOP'  
    Stop temp control (pump and heating/cooling)  
TEMP_SET_POINT = 'OUT_SP_00_'  
    Define temperature set point
```

```
UPPER_TEMP = 'OUT_SP_04_'
```

Define upper temp limit

```
build_str(param: str = "", terminator: str = '\r\n')
```

Build a command string for sending to the device

Parameters

- **param** – Command's parameter given as string
- **terminator** – Command's terminator

Returns

Command's string with a parameter and terminator

```
exception LaudaProRp245eCommandError
```

Bases: *DeviceError*

Error raised when an error is returned upon a command.

```
class LaudaProRp245eConfig(temp_set_point_init: int | float = 20.0, pump_init: int = 6, upper_temp: int | float = 80.0, lower_temp: int | float = -55.0, com_time_out: int | float = 0, max_pump_level: int = 8, max_pr_number: int = 5, operation_mode: int | OperationModeEnum = OperationModeEnum.AUTO, control_mode: int | ExtControlModeEnum = ExtControlModeEnum.INTERNAL)
```

Bases: *object*

Configuration for the Lauda RP245E circulation chiller.

```
class ExtControlModeEnum(value)
```

Bases: *IntEnum*

Source for definition of external, controlled temperature (option 2, 3 and 4 are not available with current configuration of the Lauda RP245E, add-on hardware would required)

ANALOG = 2

ETH = 5

EXPT100 = 1

INTERNAL = 0

SERIAL = 3

USB = 4

```
class OperationModeEnum(value)
```

Bases: *IntEnum*

Operation Mode (Cooling OFF/Cooling On/AUTO - set to AUTO)

AUTO = 2

Automatically select heating/cooling

COOLOFF = 0

COOLON = 1

clean_values() → *None*

com_time_out: int | float = 0

Communication time out (0 = OFF)

control_mode: int | ExtControlModeEnum = 0

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

lower_temp: int | float = -55.0

Lower temperature limit (safe for Galden HT135 cooling liquid)

max_pr_number: int = 5

Maximum number of ramp programs that can be stored in the memory of the chiller

max_pump_level: int = 8

Highest pump level of the chiller

operation_mode: int | OperationModeEnum = 2

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

pump_init: int = 6

Default pump Level

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

temp_set_point_init: int | float = 20.0

Default temperature set point

upper_temp: int | float = 80.0

Upper temperature limit (safe for Galden HT135 cooling liquid)

class LaudaProRp245eTcpCommunication(*configuration*)Bases: *Tcp*

Implements the Communication Protocol for Lauda PRO RP245E TCP connection.

close() → *None*

Close the Lauda PRO RP245E TCP connection.

static config_cls() → type[*LaudaProRp245eTcpCommunicationConfig*]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

open() → *None*

Open the Lauda PRO RP245E TCP connection.

Raises*LaudaProRp245eCommandError* – if the connection fails.**query_command(*command*: LaudaProRp245eCommand, *param*: str = '') → str**Send and receive function. E.g. to be used when setting/changing device setting. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None**read() → str**

Receive value function. :return: reply from device as a string, the terminator, as well as the 'OK' stripped from the reply to make it directly useful as a value (e.g. in case the internal bath temperature is requested)

write_command(*command*: LaudaProRp245eCommand, *param*: str = '') → *None*Send command function. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None**class LaudaProRp245eTcpCommunicationConfig(*host*: str | IPv4Address | IPv6Address, *port*: int = 54321, *bufsize*: int = 1024, *wait_sec_pre_read_or_write*: int | float = 0.005, *terminator*: str = '\r\n')**Bases: *TcpCommunicationConfig*Configuration dataclass for *LaudaProRp245eTcpCommunication*.**clean_values() → *None*****force_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

terminator: str = '\r\n'

The terminator character

wait_sec_pre_read_or_write: int | float = 0.005

Delay time between commands in seconds


Module contents

Device class for controlling a Lauda PRO RP245E, circulation chiller over TCP.

hvl_ccb.dev.mbw973

Submodules

hvl_ccb.dev.mbw973.mbw973

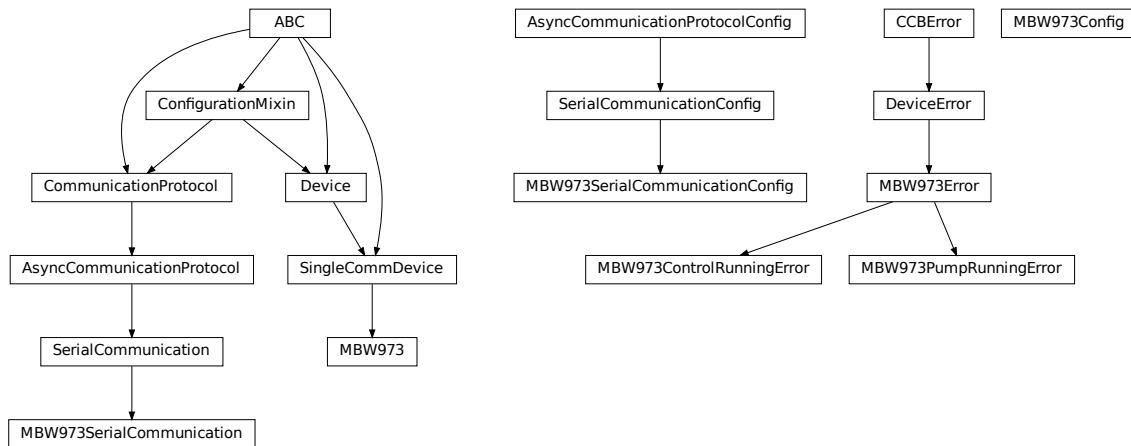

Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

```
class MBW973(com, dev_config=None)
```

Bases: *SingleCommDevice*

MBW 973 dew point mirror device class.

```
static config_cls()
```

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

```
is_done() → bool
```

Poll status of the dew point mirror and return True, if all measurements are done.

Returns

True, if all measurements are done; False otherwise.

Raises

SerialCommunicationIOError – when communication port is not opened

```
read(cast_type: type = <class 'str'>)
```

Read value from *self.com* and cast to *cast_type*. Raises *ValueError* if read text (*str*) is not convertible to *cast_type*, e.g. to *float* or to *int*.

Returns

Read value of *cast_type* type.

```
read_float() → float
```

Convenience wrapper for *self.read()*, with typing hint for return value.

Returns

Read *float* value.

```
read_int() → int
```

Convenience wrapper for *self.read()*, with typing hint for return value.

Returns

Read *int* value.

```
read_measurements() → dict[str, float]
```

Read out measurement values and return them as a dictionary.

Returns

Dictionary with values.

Raises

SerialCommunicationIOError – when communication port is not opened

```
set_measuring_options(humidity: bool = True, sf6_purity: bool = False) → None
```

Send measuring options to the dew point mirror.

Parameters

- **humidity** – Perform humidity test or not?
- **sf6_purity** – Perform SF6 purity test or not?

Raises

SerialCommunicationIOError – when communication port is not opened

start() → *None*

Start this device. Opens the communication protocol and retrieves the set measurement options from the device.

Raises

SerialCommunicationIOError – when communication port cannot be opened.

start_control() → *None*

Start dew point control to acquire a new value set.

Raises

SerialCommunicationIOError – when communication port is not opened

stop() → *None*

Stop the device. Closes also the communication protocol.

write(*value*) → *None*

Send *value* to *self.com*.

Parameters

value – Value to send, converted to *str*.

Raises

SerialCommunicationIOError – when communication port is not opened

class MBW973Config(*polling_interval: int | float = 2*)

Bases: *object*

Device configuration dataclass for MBW973.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

```
polling_interval: int | float = 2
```

Polling period for `is_done` status queries [in seconds].

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
exception MBW973ControlRunningError
```

Bases: `MBW973Error`

Error indicating there is still a measurement running, and a new one cannot be started.

```
exception MBW973Error
```

Bases: `DeviceError`

General error with the MBW973 dew point mirror device.

```
exception MBW973PumpRunningError
```

Bases: `MBW973Error`

Error indicating the pump of the dew point mirror is still recovering gas, unable to start a new measurement.

```
class MBW973SerialCommunication(configuration)
```

Bases: `SerialCommunication`

Specific communication protocol implementation for the MBW973 dew point mirror. Already predefines device-specific protocol parameters in config.

```
static config_cls()
```

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class MBW973SerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, port:
                                         Optional[str] = None, baudrate: int = 9600, parity: Union[str,
                                         hvl_ccb.comm.serial.SerialCommunicationParity] =
                                         <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                         <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                         <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                         Union[int, float] = 3)
```

Bases: `SerialCommunicationConfig`

```
baudrate: int = 9600
```

Baudrate for MBW973 is 9600 baud

```
bytesize: int | SerialCommunicationBytesize = 8
```

One byte is eight bits long

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

MBW973 does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

stopbits: int | SerialCommunicationStopbits = 1

MBW973 does use one stop bit

terminator: bytes = b'\r'

The terminator is only CR

timeout: int | float = 3

use 3 seconds timeout as default

Module contents


Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

hvl_ccb.dev.newport

Submodules

hvl_ccb.dev.newport.newport

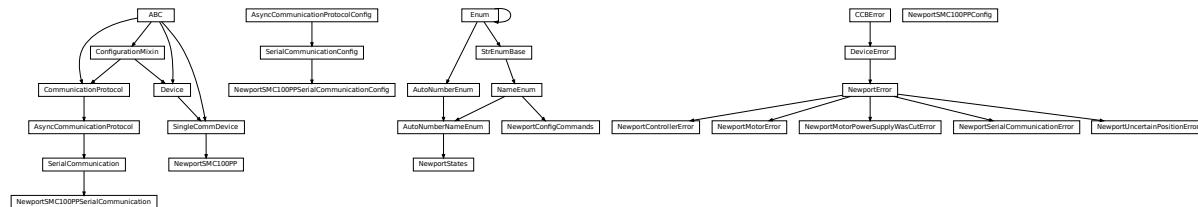

Device class for Newport SMC100PP stepper motor controller with serial communication.

The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/f/sm100-single-axis-dc-or-stepper-motion-controller>

```
class NewportConfigCommands(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *NameEnum*

Commands predefined by the communication protocol of the SMC100PP

```
AC = 'acceleration'
BA = 'backlash_compensation'
BH = 'hysteresis_compensation'
FRM = 'micro_step_per_full_step_factor'
FRS = 'motion_distance_per_full_step'
HT = 'home_search_type'
JR = 'jerk_time'
OH = 'home_search_velocity'
OT = 'home_search_timeout'
QIL = 'peak_output_current_limit'
SA = 'rs485_address'
SL = 'negative_software_limit'
SR = 'positive_software_limit'
VA = 'velocity'
VB = 'base_velocity'
ZX = 'stage_configuration'
```

exception NewportControllerError

Bases: *NewportError*

Error with the Newport controller.

exception NewportError

Bases: *DeviceError*

General Error for Newport Device

exception NewportMotorError

Bases: *NewportError*

Error with the Newport motor.

exception NewportMotorPowerSupplyWasCutError

Bases: *NewportError*

Error with the Newport motor after the power supply was cut and then restored, without interrupting the communication with the controller.

class NewportSMC100PP(*com, dev_config=None*)

Bases: *SingleCommDevice*

Device class of the Newport motor controller SMC100PP

class MotorErrors(*value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None*)

Bases: *Enum*

Possible motor errors reported by the motor during `get_state()`.

DC_VOLTAGE_TOO_LOW = 3

FOLLOWING_ERROR = 6

HOMING_TIMEOUT = 5

NED_END_OF_TURN = 11

OUTPUT_POWER_EXCEEDED = 2

PEAK_CURRENT_LIMIT = 9

POS_END_OF_TURN = 10

RMS_CURRENT_LIMIT = 8

SHORT_CIRCUIT = 7

WRONG_ESP_STAGE = 4

class StateMessages(*value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None*)

Bases: *Enum*

Possible messages returned by the controller on `get_state()` query.

CONFIG = '14'

DISABLE_FROM_JOGGING = '3E'

```

DISABLE_FROM_MOVING = '3D'
DISABLE_FROM_READY = '3C'
HOMING_FROM_RS232 = '1E'
HOMING_FROM_SMC = '1F'
JOGGING_FROM_DISABLE = '47'
JOGGING_FROM_READY = '46'
MOVING = '28'
NO_REF_ESP_STAGE_ERROR = '10'
NO_REF_FROM_CONFIG = '0C'
NO_REF_FROM_DISABLED = '0D'
NO_REF_FROM_HOMING = '0B'
NO_REF_FROM_JOGGING = '11'
NO_REF_FROM_MOVING = '0F'
NO_REF_FROM_READY = '0E'
NO_REF_FROM_RESET = '0A'
READY_FROM_DISABLE = '34'
READY_FROM_HOMING = '32'
READY_FROM_JOGGING = '35'
READY_FROM_MOVING = '33'

```

States

alias of [NewportStates](#)

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

exit_configuration(add: int | None = None) → None

Exit the CONFIGURATION state and go back to the NOT REFERENCED state. All configuration parameters are saved to the device's memory.

Parameters

add – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_acceleration(*add*: int | None = None) → int | float

Leave the configuration state. The configuration parameters are saved to the device's memory.

Parameters

add – controller address (1 to 31)

Returns

acceleration (preset units/s^2), value between 1e-6 and 1e12

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_controller_information(*add*: int | None = None) → str

Get information on the controller name and driver version

Parameters

add – controller address (1 to 31)

Returns

controller information

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_motor_configuration(*add*: int | None = None) → dict[str, float]

Query the motor configuration and returns it in a dictionary.

Parameters

add – controller address (1 to 31)

Returns

dictionary containing the motor's configuration

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_move_duration(*dist*: int | float, *add*: int | None = None) → float

Estimate the time necessary to move the motor of the specified distance.

Parameters

- **dist** – distance to travel
- **add** – controller address (1 to 31), defaults to self.address

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_negative_software_limit(*add: int | None = None*) → int | float

Get the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

Parameters

add – controller address (1 to 31)

Returns

negative software limit (preset units), value between -1e12 and 0

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_position(*add: int | None = None*) → float

Returns the value of the current position.

Parameters

add – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error
- **NewportUncertainPositionError** – if the position is ambiguous

get_positive_software_limit(*add: int | None = None*) → int | float

Get the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

Parameters

add – controller address (1 to 31)

Returns

positive software limit (preset units), value between 0 and 1e12

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_state(*add: int | None = None*) → *StateMessages*

Check on the motor errors and the controller state

Parameters

add – controller address (1 to 31)

Raises

- ***SerialCommunicationIOError*** – if the com is closed
- ***NewportSerialCommunicationError*** – if an unexpected answer is obtained
- ***NewportControllerError*** – if the controller reports an error
- ***NewportMotorError*** – if the motor reports an error

Returns

state message from the device (member of StateMessages)

go_home(*add*: int | None = None) → None

Move the motor to its home position.

Parameters

add – controller address (1 to 31), defaults to self.address

Raises

- ***SerialCommunicationIOError*** – if the com is closed
- ***NewportSerialCommunicationError*** – if an unexpected answer is obtained
- ***NewportControllerError*** – if the controller reports an error

go_to_configuration(*add*: int | None = None) → None

This method is executed during start(). It can also be executed after a reset(). The controller is put in CONFIG state, where configuration parameters can be changed.

Parameters

add – controller address (1 to 31)

Raises

- ***SerialCommunicationIOError*** – if the com is closed
- ***NewportSerialCommunicationError*** – if an unexpected answer is obtained
- ***NewportControllerError*** – if the controller reports an error

initialize(*add*: int | None = None) → None

Puts the controller from the NOT_REF state to the READY state. Sends the motor to its “home” position.

Parameters

add – controller address (1 to 31)

Raises

- ***SerialCommunicationIOError*** – if the com is closed
- ***NewportSerialCommunicationError*** – if an unexpected answer is obtained
- ***NewportControllerError*** – if the controller reports an error

move_to_absolute_position(*pos*: int | float, *add*: int | None = None) → None

Move the motor to the specified position.

Parameters

- **pos** – target absolute position (affected by the configured offset)
- **add** – controller address (1 to 31), defaults to self.address

Raises

- ***SerialCommunicationIOError*** – if the com is closed

- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

move_to_relative_position(*pos*: int | float, *add*: int | None = None) → None

Move the motor of the specified distance.

Parameters

- **pos** – distance to travel (the sign gives the direction)
- **add** – controller address (1 to 31), defaults to self.address

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

reset(*add*: int | None = None) → None

Resets the controller, equivalent to a power-up. This puts the controller back to NOT REFERENCED state, which is necessary for configuring the controller.

Parameters

- **add** – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

set_acceleration(*acc*: int | float, *add*: int | None = None) → None

Leave the configuration state. The configuration parameters are saved to the device's memory.

Parameters

- **acc** – acceleration (preset units/s^2), value between 1e-6 and 1e12
- **add** – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

set_motor_configuration(*add*: int | None = None, *config*: dict | None = None) → None

Set the motor configuration. The motor must be in CONFIG state.

Parameters

- **add** – controller address (1 to 31)
- **config** – dictionary containing the motor's configuration

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained

- *NewportControllerError* – if the controller reports an error

set_negative_software_limit(*lim: int | float, add: int | None = None*) → *None*

Set the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

Parameters

- **lim** – negative software limit (preset units), value between -1e12 and 0
- **add** – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

set_positive_software_limit(*lim: int | float, add: int | None = None*) → *None*

Set the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

Parameters

- **lim** – positive software limit (preset units), value between 0 and 1e12
- **add** – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

start()

Opens the communication protocol and applies the config.

Raises

- *SerialCommunicationIOError* – when communication port cannot be opened

stop() → *None*

Stop the device. Close the communication protocol.

stop_motion(*add: int | None = None*) → *None*

Stop a move in progress by decelerating the positioner immediately with the configured acceleration until it stops. If a controller address is provided, stops a move in progress on this controller, else stops the moves on all controllers.

Parameters

- **add** – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

wait_until_motor_initialized(*add*: int | None = None) → None

Wait until the motor leaves the HOMING state (at which point it should have arrived to the home position).

Parameters

add – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

```
class NewportSMC100PPConfig(address: int = 1, user_position_offset: int | float = 23.987, screw_scaling: int | float = 1, exit_configuration_wait_sec: int | float = 5, move_wait_sec: int | float = 1, acceleration: int | float = 10, backlash_compensation: int | float = 0, hysteresis_compensation: int | float = 0.015, micro_step_per_full_step_factor: int = 100, motion_distance_per_full_step: int | float = 0.01, home_search_type: int | HomeSearch = HomeSearch.HomeSwitch, jerk_time: int | float = 0.04, home_search_velocity: int | float = 4, home_search_timeout: int | float = 27.5, home_search_polling_interval: int | float = 1, peak_output_current_limit: int | float = 0.4, rs485_address: int = 2, negative_software_limit: int | float = -23.5, positive_software_limit: int | float = 25, velocity: int | float = 4, base_velocity: int | float = 0, stage_configuration: int | EspStageConfig = EspStageConfig.EnableEspStageCheck)
```

Bases: object

Configuration dataclass for the Newport motor controller SMC100PP.

```
class EspStageConfig(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: IntEnum

Different configurations to check or not the motor configuration upon power-up.

DisableEspStageCheck = 1

EnableEspStageCheck = 3

UpdateEspStageInfo = 2

```
class HomeSearch(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: IntEnum

Different methods for the motor to search its home position during initialization.

CurrentPosition = 1

EndOfRunSwitch = 4

EndOfRunSwitch_and_Index = 3

HomeSwitch = 2

HomeSwitch_and_Index = 0

```
acceleration: int | float = 10
```

```
address: int = 1
backlash_compensation: int | float = 0
base_velocity: int | float = 0
clean_values()
exit_configuration_wait_sec: int | float = 5
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
home_search_polling_interval: int | float = 1
```

```
home_search_timeout: int | float = 27.5
```

```
home_search_type: int | HomeSearch = 2
```

```
home_search_velocity: int | float = 4
```

```
hysteresis_compensation: int | float = 0.015
```

```
is_configdataclass = True
```

```
jerk_time: int | float = 0.04
```

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

```
micro_step_per_full_step_factor: int = 100
```

```
motion_distance_per_full_step: int | float = 0.01
```

```
property motor_config: dict[str, float]
```

Gather the configuration parameters of the motor into a dictionary.

Returns

dict containing the configuration parameters of the motor

```
move_wait_sec: int | float = 1
```

```
negative_software_limit: int | float = -23.5
```

```
classmethod optional_defaults() → dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

```
peak_output_current_limit: int | float = 0.4
positive_software_limit: int | float = 25
post_force_value(fieldname, value)
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
rs485_address: int = 2
screw_scaling: int | float = 1
stage_configuration: int | EspStageConfig = 3
user_position_offset: int | float = 23.987
velocity: int | float = 4
```

class NewportSMC100PPSSerialCommunication(configuration)

Bases: *SerialCommunication*

Specific communication protocol implementation for NewportSMC100 controller. Already predefines device-specific protocol parameters in config.

```
class ControllerErrors(value=<no_arg>, names=None, module=None, qualname=None, type=None,
start=1, boundary=None)
```

Bases: *Enum*

Possible controller errors with values as returned by the device in response to sent commands.

```
ADDR_INCORRECT = 'B'
CMD_EXEC_ERROR = 'V'
CMD_NOT_ALLOWED = 'D'
CMD_NOT_ALLOWED_CC = 'X'
CMD_NOT_ALLOWED_CONFIGURATION = 'I'
CMD_NOT_ALLOWED_DISABLE = 'J'
CMD_NOT_ALLOWED_HOMING = 'L'
CMD_NOT_ALLOWED_MOVING = 'M'
CMD_NOT_ALLOWED_NOT_REFERENCED = 'H'
CMD_NOT_ALLOWED_PP = 'W'
CMD_NOT_ALLOWED_READY = 'K'
CODE_OR_ADDR_INVALID = 'A'
COM_TIMEOUT = 'S'
```

```
DISPLACEMENT_OUT_OF_LIMIT = 'G'  
EEPROM_ACCESS_ERROR = 'U'  
ESP_STAGE_NAME_INVALID = 'F'  
HOME_STARTED = 'E'  
NO_ERROR = '@'  
PARAM_MISSING_OR_INVALID = 'C'  
POSITION_OUT_OF_LIMIT = 'N'
```

check_for_error(*add: int*) → *None*

Ask the Newport controller for the last error it recorded.

This method is called after every command or query.

Parameters

add – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

query(*add: int, cmd: str, param: int | float | str | None = None*) → *str*

Send a query to the controller, read the answer, and check for errors. The prefix add+cmd is removed from the answer.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

Returns

the answer from the device without the prefix

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

query_multiple(*add: int, cmd: str, prefixes: list[str]*) → *list[str]*

Send a query to the controller, read the answers, and check for errors. The prefixes are removed from the answers.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **prefixes** – prefixes of each line expected in the answer

Returns

list of answers from the device without prefix

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

read_text() → str

Read one line of text from the serial port, and check for presence of a null char which indicates that the motor power supply was cut and then restored. The input buffer may hold additional data afterwards, since only one line is read.

This method uses *self.access_lock* to ensure thread-safety.

Returns

String read from the serial port; ‘’ if there was nothing to read.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *NewportMotorPowerSupplyWasCutError* – if a null char is read

send_command(add: int, cmd: str, param: int | float | str | None = None) → None

Send a command to the controller, and check for errors.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

send_stop(add: int) → None

Send the general stop ST command to the controller, and check for errors.

Parameters

add – the controller address (1 to 31)

Returns

ControllerErrors reported by Newport Controller

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained

```
class NewportSMC100PPSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'ascii',
                                                encoding_error_handling: str = 'replace',
                                                wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                                default_n_attempts_read_text_nonempty: int = 10,
                                                port: Optional[str] = None, baudrate: int = 57600,
                                                parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 10)
```

Bases: *SerialCommunicationConfig*

baudrate: int = 57600

Baudrate for NewportSMC100 controller is 57600 baud

bytesize: int | SerialCommunicationBytesize = 8

NewportSMC100 controller uses 8 bits for one data byte

encoding: str = 'ascii'

use ASCII as de-/encoding, cf. the manual

encoding_error_handling: str = 'replace'

replace bytes with instead of raising utf-8 exception when decoding fails

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

NewportSMC100 controller does not use parity

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
stopbits: int | SerialCommunicationStopbits = 1
```

NewportSMC100 controller uses one stop bit

```
terminator: bytes = b'\r\n'
```

The terminator is CR/LF

```
timeout: int | float = 10
```

use 10 seconds timeout as default

```
exception NewportSerialCommunicationError
```

Bases: *NewportError*

Communication error with the Newport controller.

```
class NewportStates(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,  
boundary=None)
```

Bases: *AutoNumberNameEnum*

States of the Newport controller. Certain commands are allowed only in certain states.

CONFIG = 3

DISABLE = 6

HOMING = 2

JOGGING = 7

MOVING = 5

NO_REF = 1

READY = 4

```
exception NewportUncertainPositionError
```

Bases: *NewportError*

Error with the position of the Newport motor.

Module contents

Device class for Newport SMC100PP stepper motor controller with serial communication.


The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/f/sm100-single-axis-dc-or-stepper-motion-controller>

`hvl_ccb.dev.pfeiffer_tpg`

Submodules

`hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg`

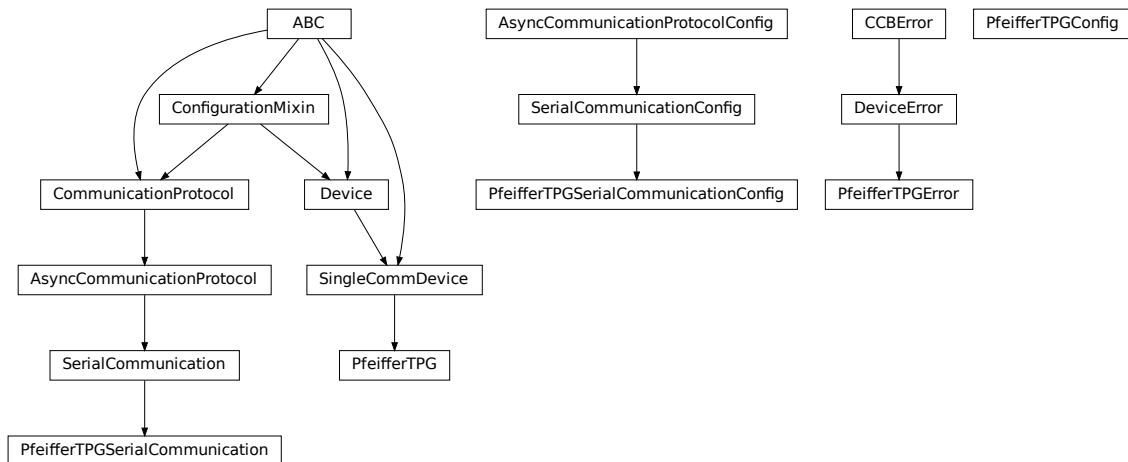

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/> measurement/activeline/controllers/

class PfeifferTPG(*com, dev_config=None*)

Bases: *SingleCommDevice*

Pfeiffer TPG control unit device class

class SensorStatus(*value*)

Bases: *IntEnum*

An enumeration.

Identification_error = 6

No_sensor = 5

Ok = 0

Overrange = 2

Sensor_error = 3

Sensor_off = 4

Underrange = 1

```
class SensorTypes(value)
```

Bases: `Enum`

An enumeration.

`CMR = 4`

`IKR = 2`

`IKR11 = 2`

`IKR9 = 2`

`IMR = 5`

`None = 7`

`PBR = 6`

`PKR = 3`

`TPR = 1`

`noSENSOR = 7`

`noSen = 7`

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

get_full_scale_mbar() → list[int | float]

Get the full scale range of the attached sensors

Returns

full scale range values in mbar, like `[0.01, 1, 0.1, 1000, 50000, 10]`

Raises

- `SerialCommunicationIOError` – when communication port is not opened
- `PfeifferTPGError` – if command fails

get_full_scale_unitless() → list[int]

Get the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

Returns

list of full scale range values, like `[0, 1, 3, 3, 2, 0]`

Raises

- `SerialCommunicationIOError` – when communication port is not opened
- `PfeifferTPGError` – if command fails

identify_sensors() → *None*

Send identification request TID to sensors on all channels.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

measure(channel: int) → tuple[str, float]

Get the status and measurement of one sensor

Parameters

channel – int channel on which the sensor is connected, with $1 \leq \text{channel} \leq \text{number_of_sensors}$

Returns

measured value as float if measurement successful, sensor status as string if not

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

measure_all() → list[tuple[str, float]]

Get the status and measurement of all sensors (this command is not available on all models)

Returns

list of measured values as float if measurements successful, and or sensor status as strings if not

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

property number_of_sensors

set_full_scale_mbar(fsr: list[int | float]) → *None*

Set the full scale range of the attached sensors (in unit mbar)

Parameters

fsr – full scale range values in mbar, for example $[0.01, 1000]$

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

set_full_scale_unitless(fsr: list[int]) → *None*

Set the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

Parameters

fsr – list of full scale range values, like $[0, 1, 3, 3, 2, 0]$

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

start() → *None*

Start this device. Opens the communication protocol, and identify the sensors.

Raises

SerialCommunicationIOError – when communication port cannot be opened

stop() → *None*

Stop the device. Closes also the communication protocol.

property unit

The pressure unit of readings is always mbar, regardless of the display unit.

class PfeifferTPGConfig(model: str | Model = Model.TPG25xA)

Bases: *object*

Device configuration dataclass for Pfeiffer TPG controllers.

class Model(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)

Bases: *NameEnum*

An enumeration.

```
TPG25xA = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}
```

```
TPGx6x = {0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8, 50000: 9}
```

is_valid_scale_reversed_str(v: str) → bool

Check if given string represents a valid reversed scale range of a model.

Parameters

v – Reversed scale range string.

Returns

True if valid, *False* otherwise.

clean_values()**force_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

```
model: str | Model = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}
```

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

exception PfeifferTPGError

Bases: *DeviceError*

Error with the Pfeiffer TPG Controller.

class PfeifferTPGSerialCommunication(configuration)

Bases: *SerialCommunication*

Specific communication protocol implementation for Pfeiffer TPG controllers. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

query(cmd: str) → str

Send a query, then read and returns the first line from the com port.

Parameters

cmd – query message to send to the device

Returns

first line read on the com

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if the device does not acknowledge the command or if the answer from the device is empty

send_command(cmd: str) → *None*

Send a command to the device and check for acknowledgement.

Parameters

cmd – command to send to the device

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if the answer from the device differs from the expected acknowledgement character ‘chr(6)’.

```
class PfeifferTPGSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Optional[str] = None, baudrate: int = 9600, parity:
                                             Union[str,
                                               hvl_ccb.comm.serial.SerialCommunicationParity] =
                                               <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                             Union[int,
                                               hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                               <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                             Union[int,
                                               hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                               <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                             Union[int, float] = 3)
```

Bases: *SerialCommunicationConfig*

baudrate: int = 9600

Baudrate for Pfeiffer TPG controllers is 9600 baud

bytesize: int | SerialCommunicationBytesize = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

Pfeiffer TPG controllers do not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
stopbits: int | SerialCommunicationStopbits = 1
Pfeiffer TPG controllers use one stop bit

terminator: bytes = b'\r\n'
The terminator is <CR><LF>

timeout: int | float = 3
use 3 seconds timeout as default
```

Module contents

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/measurement/activeline/controllers/>

hvl_ccb.dev.picotech_pt104

NOTE: PicoSDK Python wrappers already on import attempt to load the PicoSDK library; thus, the API docs can only be generated in a system with the latter installed and are by default disabled.


To build the API docs for this submodule locally edit the `docs/hvl_ccb.dev.picotech_pt104.rst` file to remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.picotech_pt104
:parts: 1

.. automodule:: hvl_ccb.dev.picotech_pt104
:members:
:undoc-members:
:show-inheritance:
```

hvl_ccb.dev.protocols

Submodules

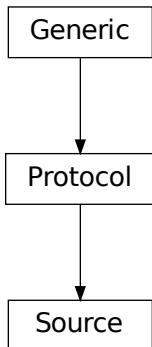hvl_ccb.dev.protocols.sources

```

class Source(*args, **kwargs)
  Bases: Protocol

  Protocol for a voltage/current source

  (Future) Implementations:
  • ea_psi9000.PSI9000
  • fug.Fug
  • heinzinger.HeinzingerPNC
  • technix.Technix

  property current: int | float
    Return the measured output current in A
  property max_current: int | float
    Maximal output current of the hardware in A, but user can reset to a lower value
  property max_voltage: int | float
    Maximal output voltage of the hardware in V, but user can reset to a lower value
  property output: bool | None
    State of the high voltage output
  property set_current: int | float
    Return the set current (may differ from actual value) in A
  property set_voltage: int | float
    Return the set voltage (may differ from actual value) in V
  property voltage: int | float
    Return the measured output voltage in V
  
```


Module contents

This module has some common protocols to interface similar device with the same methods. The benefit is to be able to switch between different devices with less effort of change the code and also to keep the same “look-and-feel”.

`hvl_ccb.dev.rs_rto1024`

Submodules

`hvl_ccb.dev.rs_rto1024.rs_rto1024`

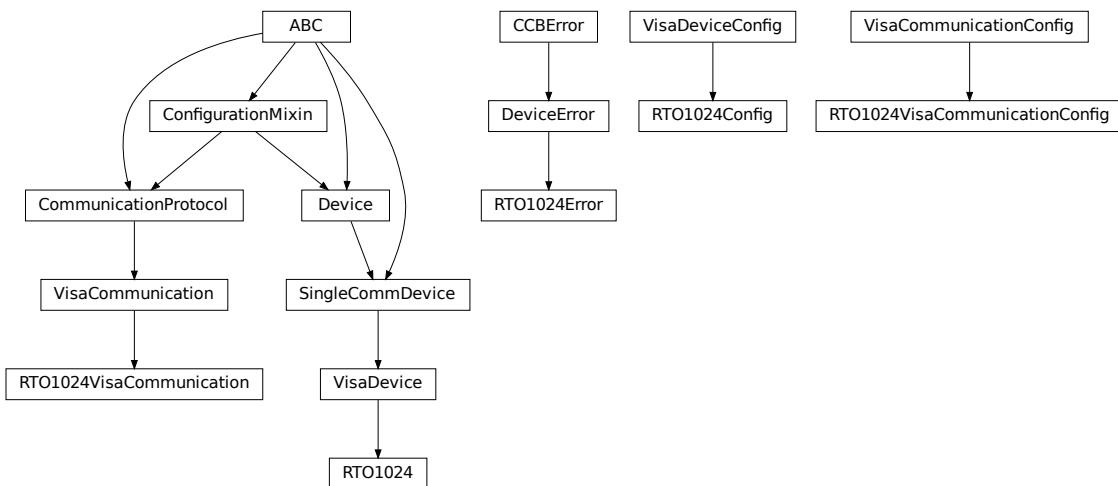

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

```
class RT01024(com: RTO1024VisaCommunication | RTO1024VisaCommunicationConfig | dict, dev_config: RTO1024Config | dict)
```

Bases: `VisaDevice`

Device class for the Rhode & Schwarz RTO 1024 oscilloscope.

```
class TriggerModes(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: `AutoNameEnum`

Enumeration for the three available trigger modes.

`AUTO = 1`

`FREERUN = 3`

`NORMAL = 2`

`classmethod names()`

Returns a list of the available trigger modes. :return: list of strings

activate_measurements(meas_n: int, source: str, measurements: list[str], category: str = 'AMPTime')

Activate the list of ‘measurements’ of the waveform ‘source’ in the measurement box number ‘meas_n’. The list ‘measurements’ starts with the main measurement and continues with additional measurements of the same ‘category’.

Parameters

- **meas_n** – measurement number 1..8
- **source** – measurement source, for example C1W1
- **measurements** – list of measurements, the first one will be the main measurement.
- **category** – the category of measurements, by default AMPTime

backup_waveform(filename: str) → None

Backup a waveform file from the standard directory specified in the device configuration to the standard backup destination specified in the device configuration. The filename has to be specified without .bin or path.

Parameters

filename – The waveform filename without extension and path

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Return the default communication protocol for this device type, which is VisaCommunication.

Returns

the VisaCommunication class

file_copy(source: str, destination: str) → None

Copy a file from one destination to another on the oscilloscope drive. If the destination file already exists, it is overwritten without notice.

Parameters

- **source** – absolute path to the source file on the DSO filesystem
- **destination** – absolute path to the destination file on the DSO filesystem

Raises

RT01024Error – if the operation did not complete

get_acquire_length() → float

Gets the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 … 500 [s]
- Increment: 1E-12 [s]

Returns

the time for one acquisition. Range: 250e-12 … 500 [s]

get_channel_offset(channel: int) → float

Gets the voltage offset of the indicated channel.

Parameters

channel – is the channel number (1..4)

Returns

channel offset voltage in V (value between -1 and 1)

get_channel_position(*channel: int*) → float

Gets the vertical position of the indicated channel.

Parameters

channel – is the channel number (1..4)

Returns

channel position in div (value between -5 and 5)

get_channel_range(*channel: int*) → float

Queries the channel range in V.

Parameters

channel – is the input channel (1..4)

Returns

channel range in V

get_channel_scale(*channel: int*) → float

Queries the channel scale in V/div.

Parameters

channel – is the input channel (1..4)

Returns

channel scale in V/div

get_channel_state(*channel: int*) → bool

Queries if the channel is active or not.

Parameters

channel – is the input channel (1..4)

Returns

True if active, else False

get_reference_point() → int

Gets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]

Returns

the reference in %

get_repetitions() → int

Get the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- *RST = 1

Returns

the number of waveforms to acquire

get_timestamps() → list[float]

Gets the timestamps of all recorded frames in the history and returns them as a list of floats.

Returns

list of timestamps in [s]

Raises

RTO1024Error – if the timestamps are invalid

list_directory(path: str) → list[tuple[str, str, int]]

List the contents of a given directory on the oscilloscope filesystem.

Parameters

path – is the path to a folder

Returns

a list of filenames in the given folder

load_configuration(filename: str) → *None*

Load current settings from a configuration file. The filename has to be specified without base directory and ‘.dfl’ extension.

Information from the manual *ReCall* calls up the instrument settings from an intermediate memory identified by the specified number. The instrument settings can be stored to this memory using the command *SAV with the associated number. It also activates the instrument settings which are stored in a file and loaded using *MMEMemory:LOAD:STATe*.

Parameters

filename – is the name of the settings file without path and extension

local_display(state: bool) → *None*

Enable or disable local display of the scope.

Parameters

state – is the desired local display state

prepare_ultra_segmentation() → *None*

Make ready for a new acquisition in ultra segmentation mode. This function does one acquisition without ultra segmentation to clear the history and prepare for a new measurement.

read_measurement(meas_n: int, name: str) → float

Parameters

- **meas_n** – measurement number 1..8
- **name** – measurement name, for example “MAX”

Returns

measured value

run_continuous_acquisition() → *None*

Start acquiring continuously.

run_single_acquisition() → *None*

Start a single or Nx acquisition.

save_configuration(*filename*: str) → *None*

Save the current oscilloscope settings to a file. The filename has to be specified without path and '.df1' extension, the file will be saved to the configured settings directory.

Information from the manual SAVe stores the current instrument settings under the specified number in an intermediate memory. The settings can be recalled using the command *RCL with the associated number. To transfer the stored instrument settings to a file, use *MMEMory:STORe:STATe* .

Parameters

filename – is the name of the settings file without path and extension

save_waveform_history(*filename*: str, *channel*: int, *waveform*: int = 1) → *None*

Save the history of one channel and one waveform to a .bin file. This function is used after an acquisition using sequence trigger mode (with or without ultra segmentation) was performed.

Parameters

- **filename** – is the name (without extension) of the file
- **channel** – is the channel number
- **waveform** – is the waveform number (typically 1)

Raises

RT01024Error – if storing waveform times out

set_acquire_length(*timerange*: float) → *None*

Defines the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]
- *RST = 0.5 [s]

Parameters

timerange – is the time for one acquisition. Range: 250e-12 ... 500 [s]

set_channel_offset(*channel*: int, *offset*: float) → *None*

Sets the voltage offset of the indicated channel.

- Range: Dependent on the channel scale and coupling [V]
- Increment: Minimum 0.001 [V], may be higher depending on the channel scale and coupling
- *RST = 0

Parameters

- **channel** – is the channel number (1..4)
- **offset** – Offset voltage. Positive values move the waveform down, negative values move it up.

set_channel_position(*channel*: int, *position*: float) → *None*

Sets the vertical position of the indicated channel as a graphical value.

- Range: -5.0 ... 5.0 [div]
- Increment: 0.02
- *RST = 0

Parameters

- **channel** – is the channel number (1..4)
- **position** – is the position. Positive values move the waveform up, negative values move it down.

set_channel_range(*channel*: int, *v_range*: float) → *None*

Sets the voltage range across the 10 vertical divisions of the diagram. Use the command alternatively instead of set_channel_scale.

- Range for range: Depends on attenuation factors and coupling. With 1:1 probe and external attenuations and 50 input coupling, the range is 10 mV to 10 V. For 1 M input coupling, it is 10 mV to 100 V. If the probe and/or external attenuation is changed, multiply the range values by the attenuation factors.
- Increment: 0.01
- *RST = 0.5

Parameters

- **channel** – is the channel number (1..4)
- **v_range** – is the vertical range [V]

set_channel_scale(*channel*: int, *scale*: float) → *None*

Sets the vertical scale for the indicated channel. The scale value is given in volts per division.

- Range for scale: depends on attenuation factor and coupling. With 1:1 probe and external attenuations and 50 input coupling, the vertical scale (input sensitivity) is 1 mV/div to 1 V/div. For 1 M input coupling, it is 1 mV/div to 10 V/div. If the probe and/or external attenuation is changed, multiply the values by the attenuation factors to get the actual scale range.
- Increment: 1e-3
- *RST = 0.05

See also: set_channel_range

Parameters

- **channel** – is the channel number (1..4)
- **scale** – is the vertical scaling [V/div]

set_channel_state(*channel*: int, *state*: bool) → *None*

Switches the channel signal on or off.

Parameters

- **channel** – is the input channel (1..4)
- **state** – is True for on, False for off

set_reference_point(*percentage*: int) → *None*

Sets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 ... 100 [%]
- Increment: 1 [%]
- *RST = 50 [%]

Parameters

percentage – is the reference in %

set_repetitions(*number: int*) → *None*

Set the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 ... 16777215
- Increment: 10
- *RST = 1

Parameters

number – is the number of waveforms to acquire

set_trigger_level(*channel: int, level: float, event_type: int = 1*) → *None*

Sets the trigger level for the specified event and source.

- Range: -10 to 10 V
- Increment: 1e-3 V
- *RST = 0 V

Parameters

- **channel** – indicates the trigger source.
 - 1..4 = channel 1 to 4, available for all event types 1..3
 - 5 = external trigger input on the rear panel for analog signals, available for A-event type = 1
 - 6..9 = not available
- **level** – is the voltage for the trigger level in [V].
- **event_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

set_trigger_mode(*mode: str | TriggerModes*) → *None*

Sets the trigger mode which determines the behavior of the instrument if no trigger occurs.

Parameters

mode – is either auto, normal, or freerun.

Raises

RTO1024Error – if an invalid triggermode is selected

set_trigger_source(*channel: int, event_type: int = 1*) → *None*

Set the trigger (Event A) source channel.

Parameters

- **channel** – is the channel number (1..4)
- **event_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

start() → *None*

Start the RTO1024 oscilloscope and bring it into a defined state and remote mode.

stop() → *None*

Stop the RTO1024 oscilloscope, reset events and close communication. Brings back the device to a state where local operation is possible.

stop_acquisition() → *None*

Stop any acquisition.

```
class RT01024Config(waveforms_path: str, settings_path: str, backup_path: str, spoll_interval: int | float = 0.5,
                      spoll_start_delay: int | float = 2, command_timeout_seconds: int | float = 60,
                      wait_sec_short_pause: int | float = 0.1, wait_sec_enable_history: int | float = 1,
                      wait_sec_post_acquisition_start: int | float = 2)
```

Bases: *VisaDeviceConfig*, *_RT01024ConfigDefaultsBase*, *_RT01024ConfigBase*

Configdataclass for the RTO1024 device.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

exception RT01024Error

Bases: *DeviceError*

class RT01024VisaCommunication(*configuration*)

Bases: *VisaCommunication*

Specialization of VisaCommunication for the RTO1024 oscilloscope

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class RT01024VisaCommunicationConfig(host: str | IPv4Address | IPv6Address, interface_type: str | InterfaceType = InterfaceType.TCPIP_INSTR, board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = '')
```

Bases: *VisaCommunicationConfig*

Configuration dataclass for VisaCommunication with specifications for the RTO1024 device class.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

interface_type: str | *InterfaceType* = 2

Interface type of the VISA connection, being one of *InterfaceType*.

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

Module contents

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

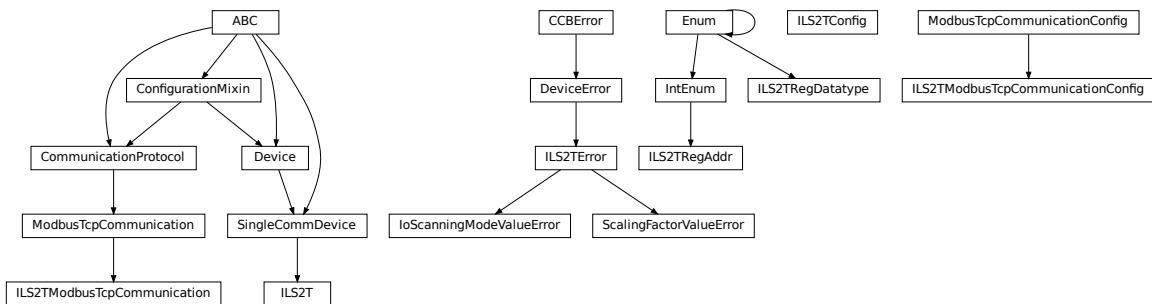hvl_ccb.dev.se_il2s**Submodules****hvl_ccb.dev.se_il2s2t.se_il2s2t**

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

class ILS2T(*com, dev_config=None*)

Bases: *SingleCommDevice*

Schneider Electric ILS2T stepper drive class.

ACTION_JOG_VALUE = 0

The single action value for *ILS2T.Mode.JOG*

class ActionsPtp(*value*)

Bases: *IntEnum*

Allowed actions in the point to point mode (*ILS2T.Mode.PTP*).

ABSOLUTE_POSITION = 0

RELATIVE_POSITION_MOTOR = 2

RELATIVE_POSITION_TARGET = 1

**DEFAULT_IO_SCANNING_CONTROL_VALUES = {
 'action': 2, 'continue_after_stop_cu': 0,
 'disable_driver_di': 0, 'enable_driver_en': 0, 'execute_stop_sh': 0,
 'fault_reset_fr': 0, 'mode': 3, 'quick_stop_qs': 0, 'ref_16': 1500, 'ref_32':
 0, 'reset_stop_ch': 0}**

Default IO Scanning control mode values

class Mode(*value*)

Bases: *IntEnum*

ILS2T device modes

JOG = 1

PTP = 3

```
class Ref16Jog(value)
```

Bases: Flag

Allowed values for ILS2T ref_16 register (the shown values are the integer representation of the bits), all in Jog mode = 1

```
FAST = 4
```

```
NEG = 2
```

```
NEG_FAST = 6
```

```
NONE = 0
```

```
POS = 1
```

```
POS_FAST = 5
```

RegAddr

Modbus Register Adresses

alias of [ILS2TRegAddr](#)

RegDatatype

Modbus Register Datatypes

alias of [ILS2TRegDatatype](#)

```
class State(value)
```

Bases: IntEnum

State machine status values

```
ON = 6
```

```
QUICKSTOP = 7
```

```
READY = 4
```

```
static config_cls()
```

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

```
disable(log_warn: bool = True, wait_sec_max: int | None = None) → bool
```

Disable the driver of the stepper motor and enable the brake.

Note: the driver cannot be disabled if the motor is still running.

Parameters

- **log_warn** – if log a warning in case the motor cannot be disabled.
- **wait_sec_max** – maximal wait time for the motor to stop running and to disable it; by default, with *None*, use a config value

Returns

True if disable request could and was sent, *False* otherwise.

do_ioscanning_write(kwargs: int) → None**

Perform a write operation using IO Scanning mode.

Parameters

kwargs – Keyword-argument list with options to send, remaining are taken from the defaults.

enable() → None

Enable the driver of the stepper motor and disable the brake.

execute_absolute_position(position: int) → bool

Execute a absolute position change, i.e. enable motor, perform absolute position change, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

Parameters

position – absolute position of motor in user defined steps.

Returns

True if actual position is as expected, *False* otherwise.

execute_relative_step(steps: int) → bool

Execute a relative step, i.e. enable motor, perform relative steps, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

Parameters

steps – Number of steps.

Returns

True if actual position is as expected, *False* otherwise.

get_dc_volt() → float

Read the DC supply voltage of the motor.

Returns

DC input voltage.

get_error_code() → dict[int, dict[str, Any]]

Read all messages in fault memory. Will read the full error message and return the decoded values. At the end the fault memory of the motor will be deleted. In addition, reset_error is called to re-enable the motor for operation.

Returns

Dictionary with all information

get_position() → int

Read the position of the drive and store into status.

Returns

Position step value

get_status() → dict[str, int]

Perform an IO Scanning read and return the status of the motor.

Returns

dict with status information.

get_temperature() → int

Read the temperature of the motor.

Returns

Temperature in degrees Celsius.

jog_run(direction: bool = True, fast: bool = False) → *None*

Slowly turn the motor in positive direction.

jog_stop() → *None*

Stop turning the motor in Jog mode.

quickstop() → *None*

Stops the motor with high deceleration rate and falls into error state. Reset with *reset_error* to recover into normal state.

reset_error() → *None*

Resets the motor into normal state after quick stop or another error occurred.

set_jog_speed(slow: int = 60, fast: int = 180) → *None*

Set the speed for jog mode. Default values correspond to startup values of the motor.

Parameters

- **slow** – RPM for slow jog mode.
- **fast** – RPM for fast jog mode.

set_max_acceleration(rpm_minute: int) → *None*

Set the maximum acceleration of the motor.

Parameters

rpm_minute – revolution per minute per minute

set_max_deceleration(rpm_minute: int) → *None*

Set the maximum deceleration of the motor.

Parameters

rpm_minute – revolution per minute per minute

set_max_rpm(rpm: int) → *None*

Set the maximum RPM.

Parameters

rpm – revolution per minute (0 < rpm <= RPM_MAX)

Raises

ILS2TError – if RPM is out of range

set_ramp_type(ramp_type: int = -1) → *None*

Set the ramp type. There are two options available:

0: linear ramp -1: motor optimized ramp

Parameters

ramp_type – 0: linear ramp | -1: motor optimized ramp

start() → *None*

Start this device.

stop() → *None*

Stop this device. Disables the motor (applies brake), disables access and closes the communication protocol.

user_steps(steps: int = 16384, revolutions: int = 1) → *None*

Define steps per revolution. Default is 16384 steps per revolution. Maximum precision is 32768 steps per revolution.

Parameters

- **steps** – number of steps in *revolutions*.
- **revolutions** – number of revolutions corresponding to *steps*.

write_absolute_position(position: int) → *None*

Write instruction to turn the motor until it reaches the absolute position. This function does not enable or disable the motor automatically.

Parameters

position – absolute position of motor in user defined steps.

write_relative_step(steps: int) → *None*

Write instruction to turn the motor the relative amount of steps. This function does not enable or disable the motor automatically.

Parameters

steps – Number of steps to turn the motor.

```
class ILS2TConfig(rpm_max_init: Integral = 1500, wait_sec_post_enable: int | float = 1,
                    wait_sec_max_disable: int | float = 10, wait_sec_post_cannot_disable: int | float = 1,
                    wait_sec_post_relative_step: int | float = 2, wait_sec_post_absolute_position: int | float = 2)
```

Bases: `object`

Configuration for the ILS2T stepper motor device.

clean_values()**force_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

rpm_max_init: Integral = 1500

initial maximum RPM for the motor, can be set up to 3000 RPM. The user is allowed to set a new max RPM at runtime using `ILS2T.set_max_rpm()`, but the value must never exceed this configuration setting.

wait_sec_max_disable: int | float = 10

wait_sec_post_absolute_position: int | float = 2

wait_sec_post_CANNOT_DISABLE: int | float = 1

wait_sec_post_enable: int | float = 1

wait_sec_post_relative_step: int | float = 2

exception ILS2TError

Bases: `DeviceError`

Error to indicate problems with the SE ILS2T stepper motor.

class ILS2TModbusTcpCommunication(configuration)

Bases: `ModbusTcpCommunication`

Specific implementation of Modbus/TCP for the Schneider Electric ILS2T stepper motor.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

class ILS2TModbusTcpCommunicationConfig(host: str | IPv4Address | IPv6Address, unit: int = 255, port: int = 502)

Bases: `ModbusTcpCommunicationConfig`

Configuration dataclass for Modbus/TCP communication specific for the Schneider Electric ILS2T stepper motor.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

unit: int = 255

The unit has to be 255 such that IO scanning mode works.

class ILS2TRegAddr(value)

Bases: IntEnum

Modbus Register Adresses for Schneider Electric ILS2T stepper drive.

ACCESS_ENABLE = 282

FLT_INFO = 15362

FLT_MEM_DEL = 15112

FLT_MEM_RESET = 15114

IO_SCANNING = 6922

JOGN_FAST = 10506

JOGN_SLOW = 10504

POSITION = 7706

RAMP_ACC = 1556

RAMP_DECEL = 1558

RAMP_N_MAX = 1554

RAMP_TYPE = 1574

SCALE = 1550

TEMP = 7200

VOLT = 7198

```
class ILS2TRegDatatype(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: `Enum`

Modbus Register Datatypes for Schneider Electric ILS2T stepper drive.

From the manual of the drive:

datatype	byte	min	max
INT8	1 Byte	-128	127
UINT8	1 Byte	0	255
INT16	2 Byte	-32_768	32_767
UINT16	2 Byte	0	65_535
INT32	4 Byte	-2_147_483_648	2_147_483_647
UINT32	4 Byte	0	4_294_967_295
BITS	just 32bits	N/A	N/A

`INT32 = (-2147483648, 2147483647)`

`is_in_range(value: int) → bool`

`exception IoScanningModeValueError`

Bases: `ILS2TError`

Error to indicate that the selected IO scanning mode is invalid.

`exception ScalingFactorValueError`

Bases: `ILS2TError`

Error to indicate that a scaling factor value is invalid.


Module contents

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

`hvl_ccb.dev.sst_luminox`

Submodules

hvl_ccb.dev.sst_luminox.sst_luminox

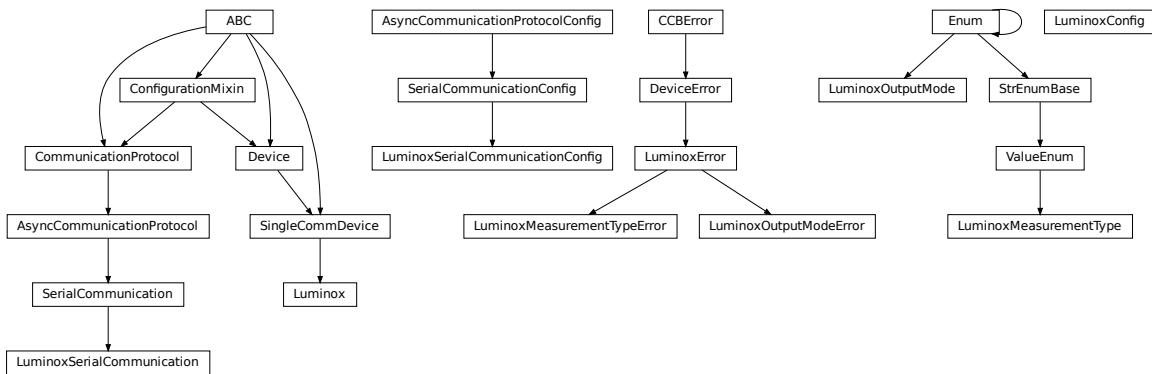

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.

Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

class Luminox(com, dev_config=None)

Bases: *SingleCommDevice*

Luminox oxygen sensor device class.

activate_output(mode: LuminoxOutputMode) → None

activate the selected output mode of the Luminox Sensor. :param mode: polling or streaming

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

query_polling(measurement: str | LuminoxMeasurementType) → dict[str | LuminoxMeasurementType, float | int | str] | float | int | str

Query a value or values of Luminox measurements in the polling mode, according to a given measurement type.

Parameters

measurement – type of measurement

Returns

value of requested measurement

Raises

- **ValueError** – when a wrong key for LuminoxMeasurementType is provided
- **LuminoxOutputModeError** – when polling mode is not activated
- **LuminoxMeasurementTypeError** – when expected measurement value is not read

read_streaming() → dict[str | *LuminoxMeasurementType*, float | int | str]

Read values of Luminox in the streaming mode. Convert the single string into separate values.

Returns

dictionary with *LuminoxMeasurementType.all_measurements_types()* keys and accordingly type-parsed values.

Raises

- **LuminoxOutputModeError** – when streaming mode is not activated
- **LuminoxMeasurementTypeError** – when any of expected measurement values is not read

start() → *None*

Start this device. Opens the communication protocol.

stop() → *None*

Stop the device. Closes also the communication protocol.

class LuminoxConfig(*wait_sec_post_activate*: int | float = 0.5, *wait_sec_trials_activate*: int | float = 0.1, *nr_trials_activate*: int = 5)

Bases: object

Configuration for the SST Luminox oxygen sensor.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

nr_trials_activate: int = 5

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

```
wait_sec_post_activate: int | float = 0.5
```

```
wait_sec_trials_activate: int | float = 0.1
```

```
exception LuminoxError
```

Bases: *DeviceError*

General Error for Luminox Device.

```
class LuminoxMeasurementType(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *ValueEnum*

Measurement types for *LuminoxOutputMode.polling*.

The *all_measurements* type will read values for the actual measurement types as given in *LuminoxOutputMode.all_measurements_types()*; it parses multiple single values using regexp's for other measurement types, therefore, no regexp is defined for this measurement type.

```
all_measurements = 'A'
```

```
classmethod all_measurements_types() → tuple[LuminoxMeasurementType, ...]
```

A tuple of *LuminoxMeasurementType* enum instances which are actual measurements, i.e. not date of manufacture or software revision.

```
barometric_pressure = 'P'
```

```
property command: str
```

```
date_of_manufacture = '# 0'
```

```
parse_read_measurement_value(read_txt: str) → dict[str | LuminoxMeasurementType, float | int | str] | float | int | str
```

```
partial_pressure_o2 = '0'
```

```
percent_o2 = '%'
```

```
sensor_status = 'e'
```

```
serial_number = '# 1'
```

```
software_revision = '# 2'
```

```
temperature_sensor = 'T'
```

```
LuminoxMeasurementTypeDict
```

A typing hint for a dictionary holding *LuminoxMeasurementType* values. Keys are allowed as strings because *LuminoxMeasurementType* is of a *StrEnumBase* type.

```
exception LuminoxMeasurementTypeError
```

Bases: *LuminoxError*

Wrong measurement type for requested data

LuminoxMeasurementTypeValue

A typing hint for all possible LuminoxMeasurementType values as read in either streaming mode or in a polling mode with *LuminoxMeasurementType.all_measurements*.

Beware: has to be manually kept in sync with *LuminoxMeasurementType* instances *cast_type* attribute values.

alias of Union[float, int, str]

class LuminoxOutputMode(value)

Bases: Enum

output mode.

polling = 1

streaming = 0

exception LuminoxOutputModeError

Bases: *LuminoxError*

Wrong output mode for requested data

class LuminoxSerialCommunication(configuration)

Bases: *SerialCommunication*

Specific communication protocol implementation for the SST Luminox oxygen sensor. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

class LuminoxSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',

encoding_error_handling: str = 'strict',

wait_sec_read_text_nonempty: Union[int, float] = 0.5,

default_n_attempts_read_text_nonempty: int = 10, port:

Optional[str] = None, baudrate: int = 9600, parity: Union[str,

hvl_ccb.comm.serial.SerialCommunicationParity] =

<SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,

hvl_ccb.comm.serial.SerialCommunicationStopbits] =

<SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,

hvl_ccb.comm.serial.SerialCommunicationBytesize] =

<SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:

Union[int, float] = 3)

Bases: *SerialCommunicationConfig*

baudrate: int = 9600

Baudrate for SST Luminox is 9600 baud

bytesize: int | SerialCommunicationBytesize = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

parity: str | SerialCommunicationParity = 'N'

SST Luminox does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

stopbits: int | SerialCommunicationStopbits = 1

SST Luminox does use one stop bit

terminator: bytes = b'\r\n'

The terminator is CR LF

timeout: int | float = 3

use 3 seconds timeout as default

Module contents

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.


Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

`hvl_ccb.dev.technix`

Submodules

`hvl_ccb.dev.technix.base`

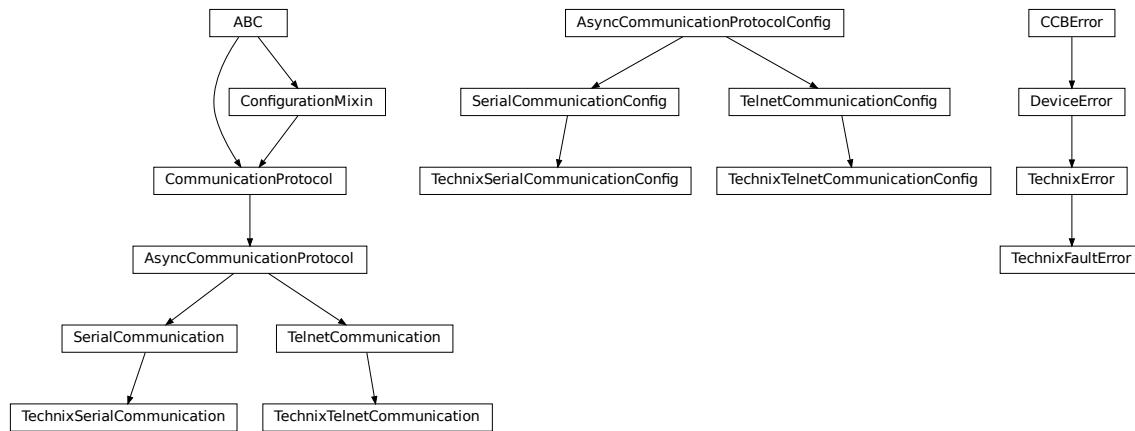

Communication and auxiliary classes for Technix

`exception TechnixError`

Bases: `DeviceError`

Technix related errors.

`exception TechnixFaultError`

Bases: `TechnixError`

Raised when the fault flag was detected while the interlock is closed

`class TechnixSerialCommunication(configuration)`

Bases: `_TechnixCommunication`, `SerialCommunication`

Serial communication for Technix

`static config_cls()`

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class TechnixSerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: int | float = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, port: str |
                                         None = None, baudrate: int = 9600, parity: str |
                                         SerialCommunicationParity =
                                         SerialCommunicationParity.NONE, stopbits: int | float |
                                         SerialCommunicationStopbits =
                                         SerialCommunicationStopbits.ONE, bytesize: int |
                                         SerialCommunicationBytesize =
                                         SerialCommunicationBytesize.EIGHTBITS, timeout: int | float =
                                         2)
```

Bases: `_TechnixCommunicationConfig, SerialCommunicationConfig`

Configuration for the serial communication for Technix

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

class TechnixTelnetCommunication(*configuration*)

Bases: `TelnetCommunication, _TechnixCommunication`

Telnet communication for Technix

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

```
class TechnixTelnetCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: int | float = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, host: str |
                                         IPv4Address | IPv6Address | None = None, port: int = 4660,
                                         timeout: int | float = 0.2)
```

Bases: `_TechnixCommunicationConfig`, `TelnetCommunicationConfig`

Configuration for the telnet communication for Technix

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

port: int = 4660


Port at which Technix is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

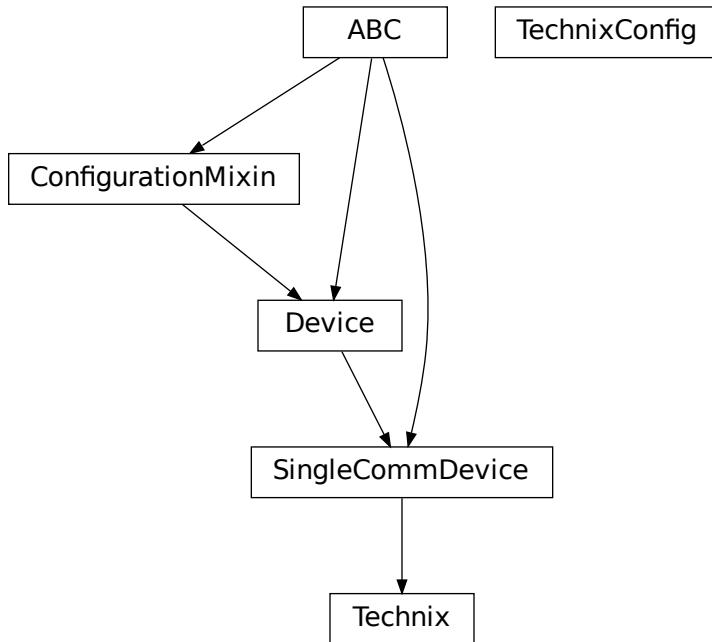hvl_ccb.dev.technix.device

its corresponding configuration class

class Technix(com, dev_config)

Bases: *SingleCommDevice*

Device class to control capacitor chargers from Technix

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property current: int | float

Actual current of the output in A

default_com_cls() → type[TechnixSerialCommunication] | type[TechnixTelnetCommunication]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

property inhibit: bool | None

Is the output of the voltage inhibited? The output stage can still be active.

property is_started: bool

Is the device started?

The device class *Technix* and

its corresponding configuration class

class Technix(com, dev_config)

Bases: *SingleCommDevice*

Device class to control capacitor chargers from Technix

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

property current: int | float

Actual current of the output in A

default_com_cls() → type[TechnixSerialCommunication] | type[TechnixTelnetCommunication]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

property inhibit: bool | None

Is the output of the voltage inhibited? The output stage can still be active.

property is_started: bool

Is the device started?

```
property max_current: int | float
    Maximal output current of the hardware in A
property max_voltage: int | float
    Maximal output voltage of the hardware in V
property open_interlock: bool | None
    Is the interlock open? (in safe mode)
property output: bool | None
    State of the high voltage output
query_status(*, _retry: bool = False)
    Query the status of the device.

Returns
    This function returns nothing

property remote: bool | None
    Is the device in remote control mode?
property set_current: int | float
    Return the set current (may differ from actual value) in A
property set_voltage: int | float
    Return the set voltage (may differ from actual value) in V
start()
    Start the device and set it into the remote controllable mode. The high voltage is turn off, and the status poller is started.
property status: _Status | None
    The status of the device with the different states as sub-fields
stop()
    Stop the device. The status poller is stopped and the high voltage output is turn off.
property voltage: int | float
    Actual voltage at the output in V
property voltage_regulation: bool | None
    Status if the output is in voltage regulation mode (or current regulation)
class TechnixConfig(communication_channel: type[hvl_ccb.dev.technix.base.TechnixSerialCommunication] |
                    type[hvl_ccb.dev.technix.base.TechnixTelnetCommunication], max_voltage: int | float,
                    max_current: int | float, polling_interval_sec: int | float = 4, post_stop_pause_sec: int | float = 1, register_pulse_time: int | float = 0.1, read_output_while_polling: bool = False)
Bases: object
clean_values()
    Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.
communication_channel: type[TechnixSerialCommunication] |
type[TechnixTelnetCommunication]
    communication channel between computer and Technix
```

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

max_current: int | float

Maximal Output current

max_voltage: int | float

Maximal Output voltage

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

polling_interval_sec: int | float = 4

Polling interval in s to maintain to watchdog of the device

post_stop_pause_sec: int | float = 1

Time to wait after stopping the device

read_output_while_polling: bool = False

Read output voltage and current within the polling event

register_pulse_time: int | float = 0.1

Time for pulsing a register

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

Module contents

Device classes for “RS 232” and “Ethernet” interfaces, which are used to control power supplies from Technix. Manufacturer homepage: <https://www.technix-hv.com>

The regulated power supplies series and capacitor chargers series from Technix are series of low and high voltage direct current power supplies as well as capacitor chargers. The class *Technix* is tested with a CCR10KV-7,5KJ via an ethernet connection as well as a CCR15-P-2500-OP via a serial connection. Check the code carefully before using it with other devices or device series


This Python package may support the following interfaces from Technix:

- Remote Interface RS232
- Ethernet Remote Interface
- Optic Fiber Remote Interface

hvl_ccb.dev.tiepie

Submodules

hvl_ccb.dev.tiepie.base

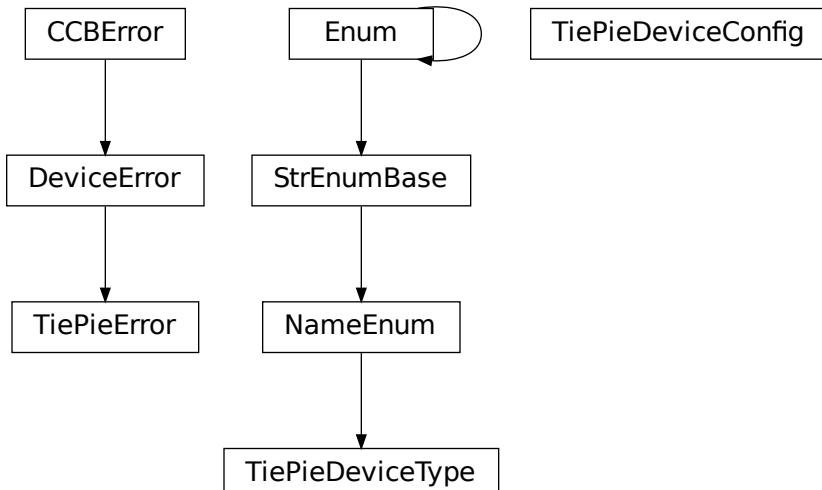

```
class TiePieDeviceConfig(serial_number: int, require_block_measurement_support: bool = True,  
                         n_max_try_get_device: int = 10, wait_sec_retry_get_device: int | float = 1.0,  
                         is_data_ready_polling_interval_sec: int | float = 0.01)
```

Bases: object

Configuration dataclass for TiePie

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

is_data_ready_polling_interval_sec: int | float = 0.01

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

n_max_try_get_device: int = 10

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

require_block_measurement_support: bool = True

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

serial_number: int

wait_sec_retry_get_device: int | float = 1.0

class TiePieDeviceType(*value=<no_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *NameEnum*

TiePie device type.

GENERATOR = 2

OSCILLOSCOPE = 1

exception TiePieError

Bases: *DeviceError*

Error of the class TiePie

```
get_device_by_serial_number(serial_number: int, device_type: str | tuple[int, _LtpDeviceReturnType],  
    n_max_try_get_device: int = 10, wait_sec_retry_get_device: float = 1.0) →  
    _LtpDeviceReturnType
```

Open and return handle of TiePie device with a given serial number

Parameters

- **serial_number** – int serial number of the device
- **device_type** – a *TiePieDeviceType* instance containing device identifier (int number) and its corresponding class, both from *libtiepie*, or a string name of such instance
- **n_max_try_get_device** – maximal number of device list updates (int number)
- **wait_sec_retry_get_device** – waiting time in seconds between retries (int number)

Returns

Instance of a *libtiepie* device class according to the specified *device_type*

Raises

- **TiePieError** – when there is no device with given serial number
- **ValueError** – when *device_type* is not an instance of *TiePieDeviceType*

```
wrap_libtiepie_exception(func: Callable) → Callable
```

Decorator wrapper for *libtiepie* methods that use *libtiepie.library.check_last_status_raise_on_error()* calls.

Parameters

func – Function or method to be wrapped


Raises

TiePieError – instead of *LibTiePieException* or one of its subtypes.

Returns

whatever *func* returns

hvl_ccb.dev.tiepie.channel

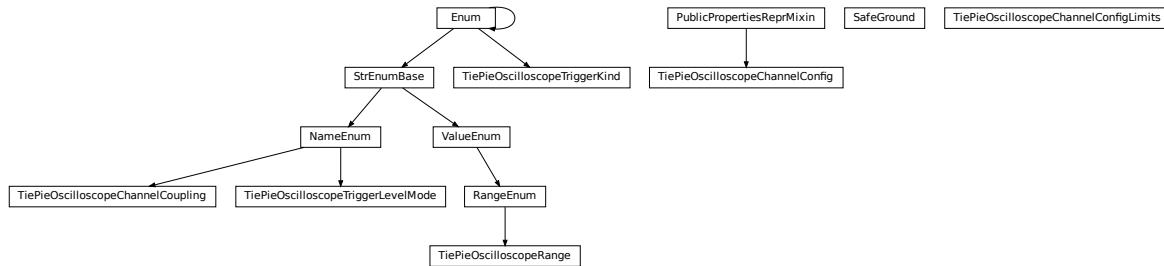

class SafeGround

Bases: object

Class that dynamically adds the `safeground_enabled` attribute getter/setter if the bound oscilloscope has the safe-ground option.

```

class TiePieOscilloscopeChannelConfig(ch_number: int, channel: OscilloscopeChannel)
Bases: PublicPropertiesReprMixin

Oscilloscope's channel configuration, with cleaning of values in properties setters as well as setting and reading them on and from the device's channel.

static clean_coupling(coupling: str | TiePieOscilloscopeChannelCoupling) →
    TiePieOscilloscopeChannelCoupling

static clean_enabled(enabled: bool) → bool

clean_input_range(input_range: float | TiePieOscilloscopeRange) → TiePieOscilloscopeRange

clean_probe_offset(probe_offset: float) → float

static clean_trigger_enabled(trigger_enabled)

clean_trigger_hysteresis(trigger_hysteresis: float) → float

static clean_trigger_kind(trigger_kind: str | TiePieOscilloscopeTriggerKind) →
    TiePieOscilloscopeTriggerKind

clean_trigger_level(trigger_level: int | float) → float

static clean_trigger_level_mode(level_mode: str | TiePieOscilloscopeTriggerLevelMode) →
    TiePieOscilloscopeTriggerLevelMode

property coupling: TiePieOscilloscopeChannelCoupling
property enabled: bool
property has_safeground: bool
    Check whether bound oscilloscope device has "safeground" option
Returns
    bool: 1=safeground available

property input_range: TiePieOscilloscopeRange
property probe_offset: float
property trigger_enabled: bool
property trigger_hysteresis: float
property trigger_kind: TiePieOscilloscopeTriggerKind
property trigger_level: float
property trigger_level_mode: TiePieOscilloscopeTriggerLevelMode

class TiePieOscilloscopeChannelConfigLimits(osc_channel: OscilloscopeChannel)
Bases: object

Default limits for oscilloscope channel parameters.

class TiePieOscilloscopeChannelCoupling(value=<no_arg>, names=None, module=None,
                                         qualname=None, type=None, start=1, boundary=None)

Bases: NameEnum

An enumeration.

```

```
ACA = 8  
ACV = 2  
DCA = 4  
DCV = 1
```

```
class TiePieOscilloscopeRange(value=<no_arg>, names=None, module=None, qualname=None,  
                               type=None, start=1, boundary=None)
```

Bases: *RangeEnum*

An enumeration.

```
EIGHTY_VOLT = 80.0
```

```
EIGHT_HUNDRED_MILLI_VOLT = 0.8
```

```
EIGHT_VOLT = 8.0
```

```
FORTY_VOLT = 40.0
```

```
FOUR_HUNDRED_MILLI_VOLT = 0.4
```

```
FOUR_VOLT = 4.0
```

```
TWENTY_VOLT = 20.0
```

```
TWO_HUNDRED_MILLI_VOLT = 0.2
```

```
TWO_VOLT = 2.0
```

```
class TiePieOscilloscopeTriggerKind(value=<no_arg>, names=None, module=None, qualname=None,  
                                      type=None, start=1, boundary=None)
```

Bases: *Enum*

An enumeration.

```
ANY = 16
```

```
FALLING = 2
```

```
RISING = 1
```

```
class TiePieOscilloscopeTriggerLevelMode(value=<no_arg>, names=None, module=None,  
                                         qualname=None, type=None, start=1, boundary=None)
```


Bases: *NameEnum*

An enumeration.

```
ABSOLUTE = 2
```

```
RELATIVE = 1
```

```
UNKNOWN = 0
```
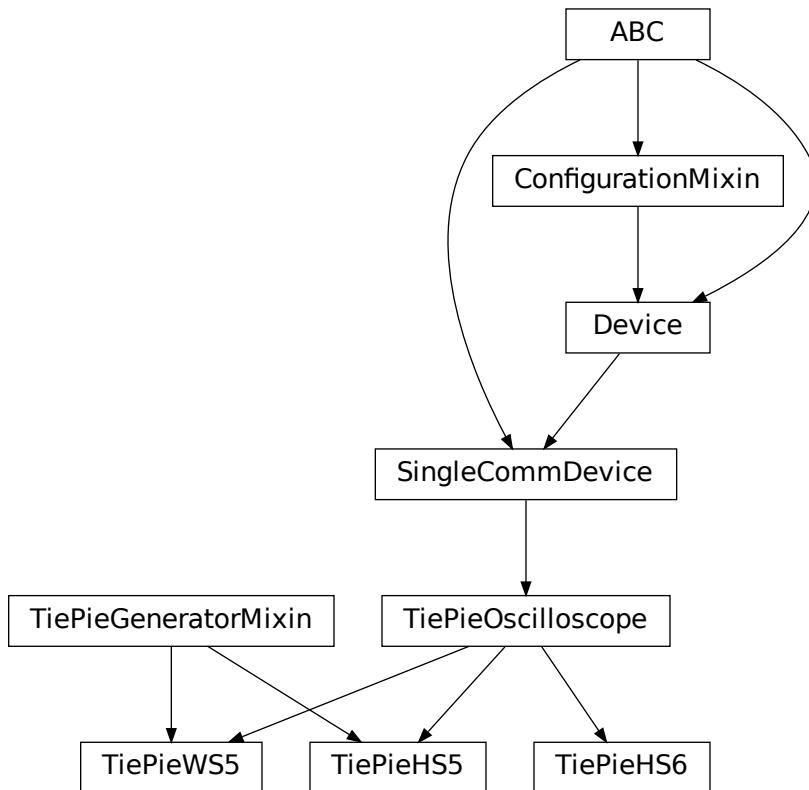
`hvl_ccb.dev.tiepie.device`

TiePie devices.

`class TiePieHS5(com, dev_config)`
 Bases: `TiePieGeneratorMixin, TiePieOscilloscope`

TiePie HS5 device.


`class TiePieHS6(com, dev_config)`
 Bases: `TiePieOscilloscope`

TiePie HS6 DIFF device.

`class TiePieWS5(com, dev_config)`
 Bases: `TiePieGeneratorMixin, TiePieOscilloscope`

TiePie WS5 device.

hvl_ccb.dev.tiepie.generator

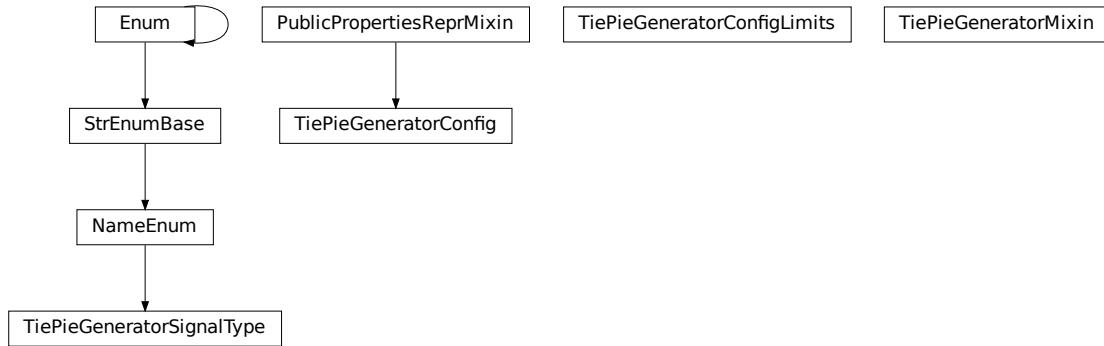

```
class TiePieGeneratorConfig(dev_gen: Generator)
    Bases: PublicPropertiesReprMixin
    Generator's configuration with cleaning of values in properties setters.

    property amplitude: float
        clean_amplitude(amplitude: float) → float
    static clean_enabled(enabled: bool) → bool
        clean_frequency(frequency: float) → float
        clean_offset(offset: float) → float
    static clean_signal_type(signal_type: int | TiePieGeneratorSignalType) → TiePieGeneratorSignalType
        clean_waveform(waveform: ndarray[Any, dtype[_ScalarType_co]]) → ndarray[Any,
            dtype[_ScalarType_co]]
    property enabled: bool
    property frequency: float
    property offset: float
    property signal_type: TiePieGeneratorSignalType
    property waveform: ndarray[Any, dtype[_ScalarType_co]] | None

class TiePieGeneratorConfigLimits(dev_gen: Generator)
    Bases: object
    Default limits for generator parameters.
```

```
class TiePieGeneratorMixin(com, dev_config)
Bases: object
TiePie Generator sub-device.
A wrapper for the libtiepie.generator.Generator class. To be mixed in with TiePieOscilloscope base class.
config_gen: TiePieGeneratorConfig | None
    Generator's dynamical configuration.

generator_start()
    Start signal generation.


generator_stop()
    Stop signal generation.

start() → None
    Start the Generator.

stop() → None
    Stop the generator.

class TiePieGeneratorSignalType(value=<no_arg>, names=None, module=None, qualname=None,
                                 type=None, start=1, boundary=None)
Bases: NameEnum
An enumeration.

ARBITRARY = 32
DC = 8
NOISE = 16
PULSE = 64
SINE = 1
SQUARE = 4
TRIANGLE = 2
UNKNOWN = 0
```

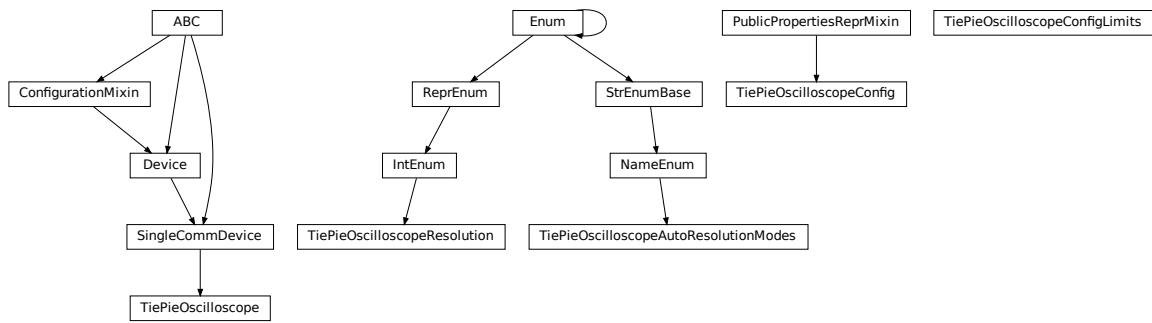hvl_ccb.dev.tiepie.oscilloscope

class TiePieOscilloscope(*com, dev_config*)

Bases: *SingleCommDevice*

TiePie oscilloscope.

A wrapper for TiePie oscilloscopes, based on the class *libtiepie.oscilloscope.Oscilloscope* with simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and its channels, including extra validation and typing hints support for configurations.

Note that, in contrast to *libtiepie* library, since all physical TiePie devices include an oscilloscope, this is the base class for all physical TiePie devices. The additional TiePie sub-devices: “Generator” and “I2CHost”, are mixed-in to this base class in subclasses.

The channels use *1..N* numbering (not *0..N-1*), as in, e.g., the Multi Channel software.

property channels_enabled: Generator[int, None, None]

Yield numbers of enabled channels.

Returns

Numbers of enabled channels

collect_measurement_data(*timeout: int | float | None = 0*) → ndarray[Any, dtype[_ScalarType_co]] | None

Try to collect the data from TiePie; return *None* if data is not ready.

Parameters

timeout – The timeout to wait until data is available. This option makes this function blocking the code. *timeout = None* blocks the code infinitely till data will be available. Per default, the *timeout* is set to *0*: The function will not block.

Returns

Measurement data of only enabled channels and time vector in a 2D-*numpy.ndarray* with float sample data; or *None* if there is no data available.

static config_cls() → type[TiePieDeviceConfig]

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

config_osc: *TiePieOscilloscopeConfig* | *None*

Oscilloscope's dynamical configuration.

config_osc_channel_dict: *dict[int, TiePieOscilloscopeChannelConfig]*

Channel configuration. A *dict* mapping actual channel number, numbered *I..N*, to channel configuration. The channel info is dynamically read from the device only on the first *start()*; beforehand the *dict* is empty.

static default_com_cls() → type[*NullCommunicationProtocol*]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

force_trigger() → *None*

Forces the TiePie to trigger with a software sided trigger event.

Return None

Raises

TiePieError – when device is not started or status of underlying device gives an error

is_measurement_data_ready() → bool

Reports if TiePie has data which is ready to collect

Returns

if the data is ready to collect.

Raises

TiePieError – when device is not started or status of underlying device gives an error

is_measurement_running() → bool

Reports if TiePie measurement is running (ready for trigger)

Returns

if a TiePie measurement is running (ready for trigger)

is_triggered() → bool

Reports if TiePie has triggered. Maybe data is not yet available. One can check with the function *is_measurement_data_ready()*.

Returns

if a trigger event occurred

static list_devices() → DeviceList

List available TiePie devices.

Returns

libtiepie up to date list of devices

property n_channels

Number of channels in the oscilloscope.

Returns

Number of channels.

start() → *None*

Start the oscilloscope.

start_measurement() → *None*

Start a measurement using set configuration.

Raises

TiePieError – when device is not started, when measurement is already running, or when status of underlying device gives an error.

stop() → *None*

Stop the oscilloscope.

stop_measurement() → *None*

Stop a measurement that is already running.

Raises

TiePieError – when device is not started, when measurement is not running, or when status of underlying device gives an error

class TiePieOscilloscopeAutoResolutionModes(*value*=<no_arg>, *names*=*None*, *module*=*None*, *qualname*=*None*, *type*=*None*, *start*=1, *boundary*=*None*)

Bases: *NameEnum*

An enumeration.

ALL = 4

DISABLED = 1

NATIVEONLY = 2

UNKNOWN = 0

class TiePieOscilloscopeConfig(*dev_osc*: Oscilloscope)

Bases: *PublicPropertiesReprMixin*

Oscilloscope's configuration with cleaning of values in properties setters.

property auto_resolution_mode: TiePieOscilloscopeAutoResolutionModes

static clean_auto_resolution_mode(*auto_resolution_mode*: int | TiePieOscilloscopeAutoResolutionModes) → TiePieOscilloscopeAutoResolutionModes

clean_pre_sample_ratio(*pre_sample_ratio*: float) → float

clean_record_length(*record_length*: int | float) → int

static clean_resolution(*resolution*: int | TiePieOscilloscopeResolution) → TiePieOscilloscopeResolution

clean_sample_rate(*sample_rate*: float) → float

clean_trigger_timeout(*trigger_timeout*: int | float | None) → float

property pre_sample_ratio: float

property record_length: int

property resolution: TiePieOscilloscopeResolution

property sample_frequency

For backwards compatibility. Use *sample_rate* instead

```
property sample_rate: float
property trigger_timeout: float | None

class TiePieOscilloscopeConfigLimits(dev_osc: Oscilloscope)
Bases: object
Default limits for oscilloscope parameters.

class TiePieOscilloscopeResolution(value=<no_arg>, names=None, module=None, qualname=None,
type=None, start=1, boundary=None)
Bases: IntEnum
An enumeration.

EIGHT_BIT = 8
FOURTEEN_BIT = 14
SIXTEEN_BIT = 16
TWELVE_BIT = 12
```

hvl_ccb.dev.tiepie.utils

PublicPropertiesReprMixin

```
class PublicPropertiesReprMixin
Bases: object
General purpose utility mixin that overwrites object representation to a one analogous to dataclass instances, but using public properties and their values instead of fields.
```

Module contents

This module is a wrapper around LibTiePie SDK devices; see <https://www.tiepie.com/en/libtiepie-sdk> .

The device classes adds simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and oscilloscope's channels. This includes extra validation and typing hints support.

Extra installation

LibTiePie SDK library is available only on Windows and on Linux.

To use this LibTiePie SDK devices wrapper:

1. install the hvl_ccb package with a tiepie extra feature:

```
$ pip install "hvl_ccb[tiepie]"
```

this will install the Python bindings for the library.

2. install the library

- on Linux: the hvl_ccb package uses the forked version

python-libtiepie-bi in version 1.1.8 which has the binaries included. No additional installation is needed anymore. * on Windows: the additional DLL is included in Python bindings package.

Troubleshooting

On a Windows system, if you encounter an OSError like this:

```
...
    self._handle = _dlopen(self._name, mode)
OSError: [WinError 126] The specified module could not be found
```


most likely the python-libtiepie package was installed in your site-packages/ directory as a python-libtiepie-*.egg file via python setup.py install or python setup.py develop command. In such case uninstall the library and re-install it using pip:

```
$ pip uninstall python-libtiepie
$ pip install python-libtiepie
```

This should create libtiepie/ folder. Alternatively, manually move the folder libtiepie/ from inside of the .egg archive file to the containing it site-packages/ directory (PyCharm's Project tool window supports reading and extracting from .egg archives).

Submodules

`hvl_ccb.dev.base`

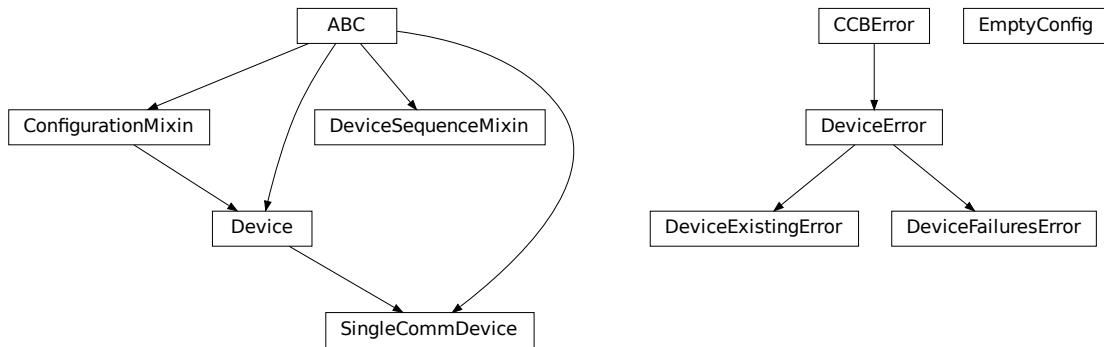

Module with base classes for devices.

class Device(`dev_config=None`)
Bases: `ConfigurationMixin`, `ABC`

Base class for devices. Implement this class for a concrete device, such as measurement equipment or voltage sources.

Specifies the methods to implement for a device.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

abstract start() → `None`

Start or restart this Device. To be implemented in the subclass.

abstract stop() → `None`

Stop this Device. To be implemented in the subclass.

exception DeviceError

Bases: `CCBError`

exception DeviceExistingError

Bases: `DeviceError`

Error to indicate that a device with that name already exists.

exception DeviceFailuresError(`failures: dict[str, Exception]`, *args)

Bases: `DeviceError`

Error to indicate that one or several devices failed.

failures: dict[str, Exception]

A dictionary of named devices failures (exceptions).

class DeviceSequenceMixin(*devices: dict[str, Device]*)

Bases: ABC

Mixin that can be used on a device or other classes to provide facilities for handling multiple devices in a sequence.

add_device(*name: str, device: Device*) → *None*

Add a new device to the device sequence.

Parameters

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

Raises

DeviceExistingError –

devices_failed_start: dict[str, Device]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices_failed_stop*.

devices_failed_stop: dict[str, Device]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices_failed_start*.

get_device(*name: str*) → *Device*

Get a device by name.

Parameters

- **name** – is the name of the device.

Returns

the device object from this sequence.

get_devices() → list[tuple[str, Device]]

Get list of name, device pairs according to current sequence.

Returns

A list of tuples with name and device each.

remove_device(*name: str*) → *Device*

Remove a device from this sequence and return the device object.

Parameters

- **name** – is the name of the device.

Returns

device object or *None* if such device was not in the sequence.

Raises

ValueError – when device with given name was not found

start() → *None*

Start all devices in this sequence in their added order.

Raises

DeviceFailuresError – if one or several devices failed to start

stop() → *None*

Stop all devices in this sequence in their reverse order.

Raises

DeviceFailuresError – if one or several devices failed to stop

class EmptyConfig

Bases: *object*

Empty configuration dataclass that is the default configuration for a Device.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

class SingleCommDevice(*com*, *dev_config=None*)

Bases: *Device*, ABC

Base class for devices with a single communication protocol.

property com

Get the communication protocol of this device.

Returns

an instance of CommunicationProtocol subtype

abstract static default_com_cls() → type[*CommunicationProtocol*]

Get the class for the default communication protocol used with this device.

Returns

the type of the standard communication protocol for this device

start() → *None*


Open the associated communication protocol.

stop() → *None*

Close the associated communication protocol.

hvl_ccb.dev.utils

hvl_ccb.dev.visa

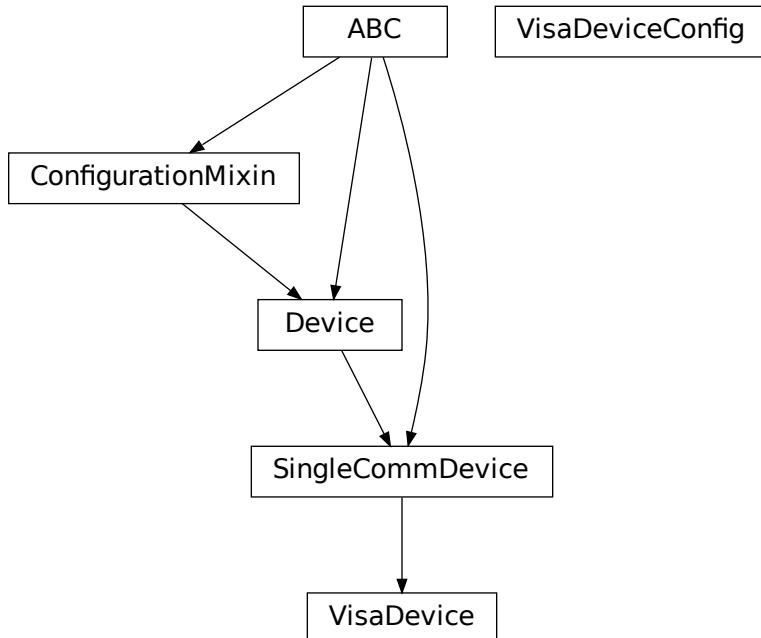

class VisaDevice(*com*: VisaCommunication | VisaCommunicationConfig | *dict*, *dev_config*: VisaDeviceConfig | *dict* | *None* = *None*)

Bases: *SingleCommDevice*

Device communicating over the VISA protocol using VisaCommunication.

static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

static default_com_cls() → type[*VisaCommunication*]

Return the default communication protocol for this device type, which is *VisaCommunication*.

Returns

the *VisaCommunication* class

get_error_queue() → str

Read out error queue and logs the error.

Returns

Error string

get_identification() → str

Queries “*IDN?” and returns the identification string of the connected device.

Returns

the identification string of the connected device

reset() → *None*

Send “*RST” and “*CLS” to the device. Typically sets a defined state.

spoll_handler()

Reads the status byte and decodes it. The status byte STB is defined in IEEE 488.2. It provides a rough overview of the instrument status.

Returns

start() → *None*

Start the *VisaDevice*. Sets up the status poller and starts it.

Returns

stop() → *None*

Stop the *VisaDevice*. Stops the polling thread and closes the communication protocol.

Returns

wait_operation_complete(timeout: float | *None* = *None*) → bool

Waits for a operation complete event. Returns after timeout [s] has expired or the operation complete event has been caught.

Parameters

timeout – Time in seconds to wait for the event; *None* for no timeout.

Returns

True, if OPC event is caught, False if timeout expired

class VisaDeviceConfig(spoll_interval: int | float = 0.5, spoll_start_delay: int | float = 2)

Bases: *_VisaDeviceConfigDefaultsBase*, *_VisaDeviceConfigBase*

Configdataclass for a VISA device.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

Module contents

Devices subpackage.

hvl_ccb.utils

Subpackages

hvl_ccb.utils.conversion

Submodules

hvl_ccb.utils.conversion.map_range

MapBitAsymRange

MapBitSymRange

MapRanges

```
class MapBitAsymRange(value: int | float, bit: int, dtype_1: ~numpy.dtype[~typing.Any] | None | type[~typing.Any] | ~numpy._typing._dtype_like._SupportsDType[~numpy.dtype[~typing.Any]] | str | tuple[~typing.Any, int] | tuple[~typing.Any, ~typing.SupportsIndex] | ~collections.abc.Sequence[~typing.SupportsIndex]] | list[~typing.Any] | ~numpy._typing._dtype_like._DTypeDict | tuple[~typing.Any, ~typing.Any] = <class 'float'>, logger=None)
```

Bases: _MapBitRange

Class to convert an asymmetric arbitrary range (0 to value) to a bit-range (0 to $2^{**\text{bit}} - 1$).

```
class MapBitSymRange(value: int | float, bit: int, dtype_1: ~numpy.dtype[~typing.Any] | None | type[~typing.Any] | ~numpy._typing._dtype_like._SupportsDType[~numpy.dtype[~typing.Any]] | str | tuple[~typing.Any, int] | tuple[~typing.Any, ~typing.SupportsIndex] | ~collections.abc.Sequence[~typing.SupportsIndex] | list[~typing.Any] | ~numpy._typing._dtype_like._DTypeDict | tuple[~typing.Any, ~typing.Any] = <class 'float'>, logger=None)
```

Bases: `_MapBitRange`

Class to convert a symmetric arbitrary range (-value to value) to a bit-range (0 to $2^{**\text{bit}} - 1$).


```
class MapRanges(range_1: tuple[int | float, int | float], range_2: tuple[int | float, int | float], dtype_1: dtype[Any] | None | type[Any] | _SupportsDType[dtype[Any]] | str | tuple[Any, int] | tuple[Any, SupportsIndex] | Sequence[SupportsIndex] | list[Any] | _DTypeDict | tuple[Any, Any], dtype_2: dtype[Any] | None | type[Any] | _SupportsDType[dtype[Any]] | str | tuple[Any, int] | tuple[Any, SupportsIndex] | Sequence[SupportsIndex] | list[Any] | _DTypeDict | tuple[Any, Any], logger=None)
```

Bases: `object`

`convert_to_range1(value, **kwargs)`

`convert_to_range2(value, **kwargs)`

[hvl_ccb.utils.conversion.sensor](#)

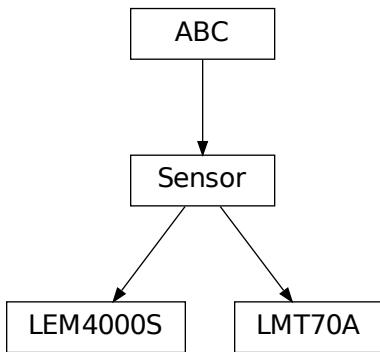

Sensors that are used by the devices implemented in the CCB

```
class LEM4000S
```

Bases: `Sensor`

`CONVERSION`

`calibration_factor`

`convert(value, **kwargs)`

`shunt`

```
class LMT70A
```

Bases: *Sensor*

Converts the output voltage (V) to the measured temperature (default °C) when using a TI Precision Analog Temperature Sensor LMT70(A)

LUT

```
convert(value, **kwargs)
```


```
temperature_unit
```

```
class Sensor
```

Bases: ABC

```
abstract convert(value, **kwargs)
```

[hvl_ccb.utils.conversion.unit](#)

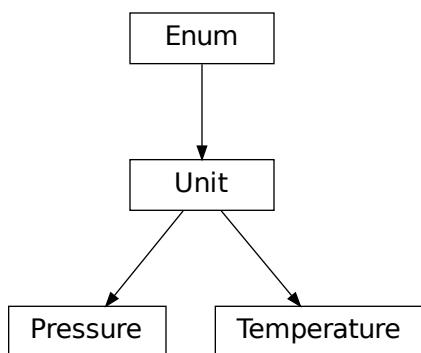

Unit conversion, within in the same group of units, for example Kelvin <-> Celsius

```
class Pressure(value)
```

Bases: *Unit*

An enumeration.

```
ATM = 'atm'
```

```
ATMOSPHERE = 'atm'
```

```
BAR = 'bar'
```

```
MILLIMETER_MERCURY = 'mmHg'
```

```
MMHG = 'mmHg'
```

```
PA = 'Pa'
```

```
PASCAL = 'Pa'
```

```

POUNDS_PER_SQUARE_INCH = 'psi'
PSI = 'psi'
TORR = 'torr'

class Temperature(value)
    Bases: Unit
    An enumeration.


    C = 'C'
    CELSIUS = 'C'
    F = 'F'
    FAHRENHEIT = 'F'
    K = 'K'
    KELVIN = 'K'

class Unit(value)
    Bases: Enum
    An enumeration.

    abstract classmethod convert(value, **kwargs)

```

[hvl_ccb.utils.conversion.utils](#)

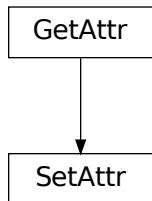

```

class GetAttr(default, name)
    Bases: object

class SetAttr(default, name, limits, absolut=False, dtype=(<class 'int'>, <class 'float'>), validator=None)
    Bases: GetAttr


convert_value_to_str(*value: ndarray[Any, dtype[_ScalarType_co]]) → list[str | list[str]]

```

Converts two sets of values to strings. This is necessary because a 0-dim array needs different treatment than a 1-dim array :param value: array of values either 0-dim or 1-dim :return:converted se

preserve_type(func)

This wrapper preserves the first order type of the input. Upto now the type of the data stored in a list, tuple, array or dict is not preserved. Integer will be converted to float!

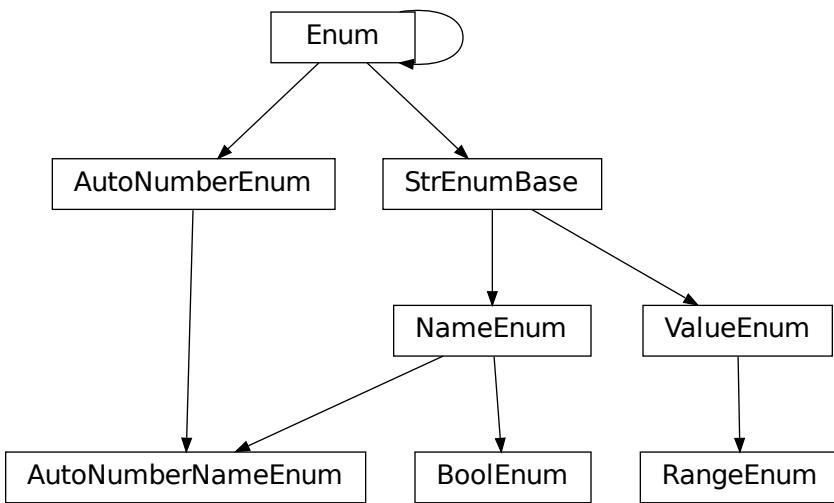Module contents**Submodules****hvl_ccb.utils.enum**

```
class AutoNumberNameEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *NameEnum*, *AutoNumberEnum*

Auto-numbered enum with names used as string representation, and with lookup and equality based on this representation.

```
class BoolEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *NameEnum*

BoolEnum inherits from NameEnum and the type of the first value is
enforced to be ‘boolean’. For bool()-operation the `__bool__` is redefined here.

```
class NameEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *StrEnumBase*

Enum with names used as string representation, and with lookup and equality based on this representation. The lookup is implemented in *StrEnumBase* with the `_missing_value_` method. The equality is also defined at this place (`__eq__`).

Use-case:

```
class E(NameEnum):
    a = 2
    b = 4

    E.a == "a"
    E.a != 2
    E.a != "2"
```

The access would be normally with E["a"], but E("a") works also. Therefore, E["a"] == E("a")

Attention: to avoid errors, best use together with *unique* enum decorator.

```
class RangeEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                boundary=None)
```

Bases: float, [ValueEnum](#)

Range enumeration inherit from ValueEnum, find suitable voltage/current/resistance input range for devices such as multimeter and oscilloscope

abstract classmethod unit() → str

Returns the Unit of the values in the enumeration. :return: the unit of the values in the enumeration in string format

```
class StrEnumBase(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
```

Bases: Enum

String representation-based equality and lookup.

```
class ValueEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                 boundary=None)
```

Bases: [StrEnumBase](#)

Enum with string representation of values used as string representation, and with lookup and equality based on this representation. Values do not need to be of type 'str', but they need to have a str-representation to enable this feature. The lookup is implemented in StrEnumBase with the `_missing_value_` method. The equality is also defined at this place (`__eq__`).

Use-case:

```
class E(ValueEnum):
    ONE = 1

    E.ONE == "1"
    E.ONE != 1
    E.ONE != "ONE"
```

The access would be normally with E(1), but E("1") works also. Therefore, E(1) == E("1")

Attention: to avoid errors, best use together with *unique* enum decorator.

hvl_ccb.utils.poller

Poller

```
class Poller(spoll_handler: Callable, polling_delay_sec: int | float = 0, polling_interval_sec: int | float = 1,  
            polling_timeout_sec: int | float | None = None)
```

Bases: `object`

Poller class wrapping `concurrent.futures.ThreadPoolExecutor` which enables passing of results and errors out of the polling thread.

is_polling() → bool

Check if device status is being polled.

Returns

True when polling thread is set and alive

start_polling() → bool

Start polling.

Returns

True if was not polling before, *False* otherwise

stop_polling() → bool

Stop polling.

Wait for until polling function returns a result as well as any exception / error that might have been raised within a thread.

Returns

True if was polling before, *False* otherwise, and last result of the polling function call.

Raises

polling function exceptions

wait_for_polling_result()

Wait for until polling function returns a result as well as any exception / error that might have been raised within a thread.

Returns

polling function result

Raises

polling function errors

hvl_ccb.utils.typing

Additional Python typing module utilities

ConvertibleTypes

Typing hint for data type that can be used in conversion

alias of Union[int, float, list[Union[int, float]], tuple[Union[int, float], ...], dict[str, Union[int, float]], ndarray[Any, dtype[_ScalarType_co]]]

Number

Typing hint auxiliary for a Python base number types: *int* or *float*.

alias of Union[int, float]

hvl_ccb.utils.validation

validate_and_resolve_host(*host*: str | IPv4Address | IPv6Address, *logger*: Logger | None = None) → str

validate_bool(*x_name*: str, *x*: object, *logger*: Logger | None = None) → None

Validate if given input *x* is a *bool*.

Parameters

- **x_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as boolean
- **logger** – logger of the calling submodule

Raises

TypeError – when the validated input does not have boolean type

validate_number(*x_name*: str, *x*: object, *limits*: tuple | None = (None, None), *number_type*: type[int | float] | tuple[type[int | float], ...] = (<class 'int'>, <class 'float'>), *logger*: ~logging.Logger | None = None) → None

Validate if given input *x* is a number of given *number_type* type, with value between given *limits[0]* and *limits[1]* (inclusive), if not *None*. For array-like objects (npt.NDArray, list, tuple, dict) it is checked if all elements are within the limits and have the correct type.

Parameters

- **x_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as number of given type within given range
- **logger** – logger of the calling submodule
- **limits** – [lower, upper] limit, with *None* denoting no limit: [-inf, +inf]
- **number_type** – expected type or tuple of types of a number, by default (*int, float*)

Raises


- **TypeError** – when the validated input does not have expected type
- **ValueError** – when the validated input has correct number type but is not within given range or has wrong input limits

validate_tcp_port(*port*: int, *logger*: Logger | None = None)

Module contents

4.1.2 Submodules

`hvl_ccb.configuration`

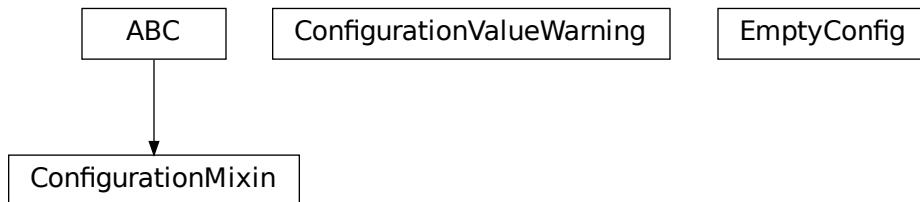

providing classes for handling configuration for communication protocols and devices.

Facilities

class ConfigurationMixin(*configuration*)

Bases: `ABC`

Mixin providing configuration to a class.

property config

ConfigDataclass property.

Returns

the configuration

abstract static config_cls()

Return the default configdataclass class.

Returns

a reference to the default configdataclass class

configuration_save_json(*path*: str) → None

Save current configuration as JSON file.

Parameters

`path` – path to the JSON file.

classmethod from_json(*filename*: str)

Instantiate communication protocol using configuration from a JSON file.

Parameters

`filename` – Path and filename to the JSON configuration

exception ConfigurationValueWarning

Bases: `UserWarning`

User warnings category for values of `@configdataclass` fields.

class EmptyConfig

Bases: `object`

Empty configuration dataclass.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

Returns

a list of strings containing all keys.

classmethod optional_defaults() → dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns

a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns

a list of strings containing all required keys.

configdataclass(direct_decoration=None, frozen=True)

Decorator to make a class a configdataclass. Types in these dataclasses are enforced. Implement a function *clean_values(self)* to do additional checking on value ranges etc.

It is possible to inherit from a configdataclass and re-decorate it with @configdataclass. In a subclass, default values can be added to existing fields. Note: adding additional non-default fields is prone to errors, since the order has to be respected through the whole chain (first non-default fields, only then default-fields).

Parameters

- **frozen** – defaults to True. False allows to later change configuration values. Attention: if configdataclass is not frozen and a value is changed, typing is not enforced anymore!

hvl_ccb.error


```
CCBError
```

Introduce a common code base error for the CCB

exception CCBError

Bases: `Exception`

hvl_ccb.experiment_manager

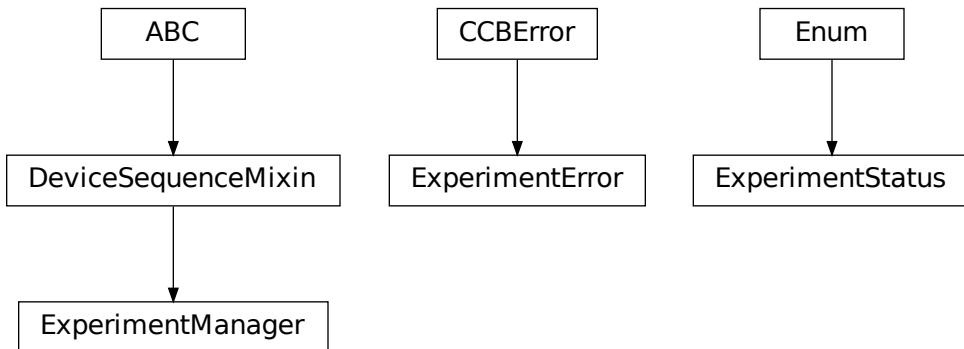

Main module containing the top level `ExperimentManager` class. Inherit from this class to implement your own experiment functionality in another project and it will help you start, stop and manage your devices.

exception ExperimentError

Bases: `CCBError`

Error to indicate that the current status of the experiment manager is on ERROR and thus no operations can be made until reset.

class ExperimentManager(*args, **kwargs)

Bases: `DeviceSequenceMixin`

Experiment Manager can start and stop communication protocols and devices. It provides methods to queue commands to devices and collect results.

add_device(name: str, device: Device) → None

Add a new device to the manager. If the experiment is running, automatically start the device. If a device with this name already exists, raise an error.

Parameters

- **name** – is the name of the device.

- **device** – is the instantiated Device object.

Raises

ExperimentError –

devices_failed_start: dict[str, Device]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices_failed_stop*.

devices_failed_stop: dict[str, Device]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices_failed_start*.

finish() → None

Stop experimental setup, stop all devices.

is_error() → bool

Returns true, if the status of the experiment manager is *error*.

Returns

True if on error, false otherwise

is_finished() → bool

Returns true, if the status of the experiment manager is *finished*.

Returns

True if finished, false otherwise

is_running() → bool

Returns true, if the status of the experiment manager is *running*.

Returns

True if running, false otherwise

run() → None

Start experimental setup, start all devices.

start() → None

Alias for ExperimentManager.run()

property status: ExperimentStatus

Get experiment status.

Returns

experiment status enum code.

stop() → None

Alias for ExperimentManager.finish()

class ExperimentStatus(value)

Bases: Enum

Enumeration for the experiment status

ERROR = 5

FINISHED = 4

FINISHING = 3

```
INITIALIZED = 0
INITIALIZING = -1
RUNNING = 2
STARTING = 1
```

4.1.3 Module contents

Top-level package for HVL Common Code Base.

CONTRIBUTING

Contributions are welcome, and they are greatly appreciated! Every little bit helps, and credit will always be given.

You can contribute in many ways:

5.1 Types of Contributions

5.1.1 Report Bugs

Report bugs at https://gitlab.com/ethz_hvl/hvl_ccb/issues.

If you are reporting a bug, please include:

- Your operating system name and version.
- Any details about your local setup that might be helpful in troubleshooting.
- Detailed steps to reproduce the bug.

5.1.2 Fix Bugs

Look through the GitLab issues for bugs. Anything tagged with “bug” and “help wanted” is open to whoever wants to implement it.

5.1.3 Implement Features

Look through the GitLab issues for features. Anything tagged with “enhancement” and “help wanted” is open to whoever wants to implement it.

5.1.4 Write Documentation

HVL Common Code Base could always use more documentation, whether as part of the official HVL Common Code Base docs, in docstrings, or even on the web in blog posts, articles, and such.

5.1.5 Submit Feedback

The best way to send feedback is to file an issue at https://gitlab.com/ethz_hvl/hvl_ccb/issues.

If you are proposing a feature:

- Explain in detail how it would work.
- Keep the scope as narrow as possible, to make it easier to implement.
- Remember that this is a volunteer-driven project, and that contributions are welcome :)

5.2 Get Started!

Ready to contribute? Here's how to set up *hvl_ccb* for local development.

1. Clone *hvl_ccb* repo from GitLab:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

2. Go into the cloned folder. Then install your virtual environment and activate it:

```
$ cd hvl_ccb
$ python -m venv .venv
$ . .venv/Scripts/activate # <-- for Windows
$ . .venv_22/bin/activate # <-- for Linux
```

3. Install the HVL-CommonCodeBase with its dependencies as well as the dependencies for development:

```
$ pip install -e .[all]
$ pip install -r requirements_dev.txt
```

4. Furthermore, it is recommended to install the git hook script shipped within the repository:

```
$ pre-commit install
```

5. After creating an Issue and Merge Request on GitLab, you can switch to your created development branch:

```
$ git switch name-of-your-bugfix-or-feature
```

Now you can make your changes locally.

6. When you're done making changes, check that your changes pass flake8, mypy, black, isort and the tests, including testing other Python versions with tox:

```
$ isort .
$ black --preview hvl_ccb/ tests/ examples/
$ flake8 hvl_ccb tests
$ mypy --show-error-codes hvl_ccb
$ python setup.py test or py.test
$ tox
```

You can also use the provided make-like shell script to run flake8 and tests:

```
$ ./make.sh black
$ ./make.sh isort
$ ./make.sh style
$ ./make.sh type
$ ./make.sh test
```

- As we want to maintain a high quality of coding style we use *black* and *isort*. This style is checked with the pipelines on gitlab.com. Ensure that your commits include only properly formatted code. One way to comply is to install and use *pre-commit*. This package includes the necessary configuration.
- Commit your changes and push your branch to GitLab:

```
$ git add .
$ git commit -m "Your detailed description of your changes."
$ git push
```

- Request a review of your merge request through the GitLab website.

5.3 Merge Request Guidelines

Before you submit a merge request, check that it meets these guidelines:

- The merge request should include tests.
- If the merge request adds functionality, the docs should be updated. Put your new functionality into a function with a docstring, and add the feature to the list in README.rst.
- The merge request should work for Python 3.9 to 3.10. Check https://gitlab.com/ethz_hvl/hvl_ccb/merge_requests and make sure that the tests pass for all supported Python versions.

5.4 Tips

- To run tests from a single file:

```
$ py.test tests/test_hvl_ccb.py
```

or a single test function:

```
$ py.test tests/test_hvl_ccb.py::test_command_line_interface
```

- If your tests are slow, profile them using the pytest-profiling plugin:

```
$ py.test tests/test_hvl_ccb.py --profile
```

or for a graphical overview (you need a SVG image viewer):

```
$ py.test tests/test_hvl_ccb.py --profile-svg
$ open prof/combined.svg
```

- To add dependency, edit appropriate *requirements variable in the setup.py file and re-run:

```
$ python setup.py develop
```

- To generate a PDF version of the Sphinx documentation instead of HTML use:

```
$ rm -rf docs/hvl_ccb.rst docs/modules.rst docs/_build && sphinx-apidoc -o docs/hvl_ccb && python -msphinx -M latexpdf docs/ docs/_build
```

This command can also be run through the make-like shell script:

```
$ ./make.sh docs-pdf
```

This requires a local installation of a LaTeX distribution, e.g. MikTeX.

5.5 Deploying

A reminder for the maintainers on how to deploy.

Make sure all your changes are committed and that all relevant MR are merged. Then switch to `devel`, update it and create `release-N.M.K` branch:

```
$ git switch devel  
$ git pull  
$ git checkout -b release-N.M.K
```

- Update copyright information (if necessary) in `docs/conf.py` and `README.rst`
- Update or create entry in `HISTORY.rst` (commit message: Update HISTORY.rst: release N.M.K).
- Update, if applicable, `AUTHORS.rst` (commit message: Update AUTHORS.rst: release N.M.K)
- Update features tables in `README.rst` file (commit message: Update README.rst: release N.M.K)
- Update API docs (commit message: Update API-docs: release N.M.K)

```
$ ./make.sh docs # windows  
$ make docs # unix-based-os
```

Commit all of the above, except for

- `docs/hvl_ccb.dev.picotech_pt104.rst`.

Before you continue revert the changes in this file.

Then run:

```
$ bumpver update --patch # possible: major / minor / patch  
$ git push --set-upstream origin release-N.M.K  
$ git push --tags
```

Go to <https://readthedocs.org/projects/hvlccb/builds/> and check if RTD docs build for the pushed tag passed.

Wait for the CI pipeline to finish successfully.

The two following commands are best executed in a WSL or Unix based OS. Run a release check:

```
$ make release-check
```

Finally, prepare and push a release:

```
$ make release
```

Merge the release branch into master and devel branches with --no-ff flag and delete the release branch:

```
$ git switch master
$ git pull
$ git merge --no-ff release-N.M.K
$ git push
$ git switch devel
$ git merge --no-ff release-N.M.K
$ git push
$ git push --delete origin release-N.M.K
$ git branch --delete release-N.M.K
```

After this you can/should clean your folder (with WSL/Unix command):

```
$ make clean
```

Finally, prepare GitLab release and cleanup the corresponding milestone:

1. go to https://gitlab.com/ethz_hvl/hvl_ccb/-/tags/, select the latest release tag, press “Edit release notes” and add the release notes (copy a corresponding entry from `HISTORY.rst` file with formatting adjusted from ReStructuredText to Markdown); press “Save changes”;
2. go to https://gitlab.com/ethz_hvl/hvl_ccb/-/releases, select the latest release, press “Edit this release” and under “Milestones” select the corresponding milestone; press “Save changes”;
3. go to https://gitlab.com/ethz_hvl/hvl_ccb/-/milestones, make sure that it is 100% complete (otherwise, create a next patch-level milestone and assign it to the ongoing Issues and Merge Requests therein); press “Close Milestone”.

**CHAPTER
SIX**

CREDITS

6.1 Active Maintainers

- Chi-Ching Hsu <hsu@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>

6.2 Authors and Contributors

- Alise Chachereau (Author)
- Maria Del (Contributor, Maintainer)
- Joseph Engelbrecht (Author)
- Raphael Faerber (Contributor)
- David Graber (Author)
- Chi-Ching Hsu (Author, Maintainer)
- Henning Janssen (Author, Maintainer)
- Henrik Menne (Author, Maintainer)
- Luca Nembrini (Contributor)
- Mikołaj Rybiński (Author, Maintainer)
- Ruben Stadler (Contributor)
- David Taylor (Author)
- Hanut Vemulapalli (Contributor)

HISTORY

7.1 0.14.4 (2023-12-22)

- Hot-fix to be compatible with the newly released version 23.12.0 of black
- Switch to src-layout
- In `tiepie`:
 - implementation property `generator_is_running` to check, whether the generator is running
 - properties `probe_offset` and `probe_gain` are not implemented and raise `NotImplementedError`

7.2 0.14.3 (2023-11-17)

- Fix Heinzinger conversion from mA to A and fix wrong docstrings
- Hot-fix to be compatible with the newly released version 23.11.0 of black
- Fix `bumpver` tag messages
- Implementation of Technix to fulfil `protocols.Source`

7.3 0.14.2 (2023-09-07)

- Change dependency to `libtiepie` with linux binaries using forked version `python-libtiepie-bi`
- Adapt `makefile` after removing `setup.py`
- Fix commit messages and tag with `bumpver`
- Hot-fix to be compatible with the newly released version 3.5.0 of `pymodbus`

7.4 0.14.1 (2023-08-21)

- Remove `setup.cfg`, `setup.py`, `mypy.ini`, `pytest.ini`, `requirements_dev.txt` and change to `pyproject.toml`
- Replace `bump2version` with `bumpver`
- Change dependency to `libtiepie` with linux binaries
- Hot-fix to be compatible with the newly released version 6.1.0 of `flake8`

7.5 0.14.0 (2023-07-28)

- **Rework of Heinzinger high voltage source control**
 - validation of input values (e.g. `voltage`-property)
 - merge `HeinzingerPNC` and `HeinzingerDI` to `Heinzinger`
 - always return values as V for voltage and A for current
 - `RangeEnum` for `number_of_recordings`
 - fulfil `protocols.Source`
 - raise Error for getter and setter (e.g. `set_current`, `get_current`...), use property instead
- Hot-fix to be compatible with the newly released version 23.7.0 of `black`
- Hot-fix to be compatible with the newly released version 3.4.0 of `pymodbus`
- **Hot-fix to be compatible with the newly released version 1.1.6 of `libtiepie`**
 - drop support for I2C
- Remove default import from `hvl_ccb.comm` and `hvl_ccb.dev` for specific communication protocols and devices

7.6 0.13.3 (2023-03-31)

- Introduce common protocol for voltage and current sources `hvl_ccb.dev.protocols.sources`
- Update code style to `black` 23.3.0 and `isort` 5.12.0
- An `_EarthingStick` of `BaseCube` is implemented as a `_Switch`
- **Code improvements for device Heinzinger**
 - use property-based instead of getter and setter
 - `DeprecationWarning` for getter and setter (e.g. `set_current`, `get_current`...)

7.7 0.13.2 (2023-03-17)

- Hot-fix to be compatible with the newly released version 3.0.0 of typeguard

7.8 0.13.1 (2023-03-03)

- **Repository maintenance**

- add the option to manually set `n_attempts_max` and `attempt_interval_sec` in query of the `SyncCommunicationProtocol`
- fix links in description for Heinzinger digital interface and universal high voltage power supplies
- keep copyright year information only in `docs/conf.py` and `README.rst`
- remove copyright year information from the files
- fix readthedocs build failed issue
- update code style to `black 23.1.0`

7.9 0.13.0 (2023-01-27)

- **Drop support for Python 3.7 and 3.8:**

- remove version dependent implementations
- changed typing acc. to PEP 585

- Un-freeze version number of dependencies and upgrade to most recent versions

7.10 0.12.3 (2022-12-27)

- **Code improvements for device cube:**

- split alarms from constants
- split errors from constants
- split earthing_stick from constants
- split support from constants

- Update code style to `black 22.12.0`

- **Smaller change of device tiepie:**

- change hard coded trigger time out value for no time out/infinite (-1) to `ltp.const.TO_INFINITY`

7.11 0.12.2 (2022-11-29)

- Move the device modules into packages
- Bugfix in `validate_number` to check the order of the limits
- **Repository maintenance:**
 - imports are sorted with `isort`
 - some `mypy` fixing and additional typing

7.12 0.12.1 (2022-10-31)

- Fix `numpy` version requirement problem
 - for Python 3.7: 1.21.6
 - for Python 3.8 and onwards: 1.23.4

7.13 0.12.0 (2022-10-17)

- Last release for Python 3.7 and 3.8
- **Repository maintenance**
 - update Labjack LJM software installer link in the pipeline
 - fix dependencies to the fixed version
 - fix `asyncua` to 0.9.95 and `pymodbus` to 2.5.3 (newer versions break the code)
 - fix PICube checker for slope as it is always positive

7.14 0.11.1 (2022-09-15)

- **Repository maintenance**
 - fix issue with `mypy` and Python 3.10.7
 - update code style to `black` 22.8.0
 - project configurations merged into `setup.cfg`
 - fix coverage indicator

7.15 0.11.0 (2022-06-22)

- New device: Fluke 884X Bench 6.5 Digit Precision Multimeter
- RangeEnum is a new enum for e.g. measurement ranges which also finds a suitable range object
- **smaller changes of device tiepie:**
 - introduce status method `is_measurement_running()` to check if the device is armed
 - introduce `stop_measurement()` to disarm the trigger of the device
 - fix bug with docs due to change of libtiepie
- NameEnum and inherited enums can only have unique entries

7.16 0.10.3 (2022-03-21)

- fix bug in the Labjack pulse feature that occurred when the start time was set to 0s
- new conversion utility to map two ranges on each other
- update CONTRIBUTING.RST
- update makefile and make.sh
- improve the mockup telnet test server

7.17 0.10.2 (2022-02-28)

- introduction of `black` as code formatter
- increase the required version of the package `aenum`
- remove device `supercube2015` - as it is no longer used
- remove unused package `openpyxl` requirement
- fix bug in highland logging
- improve handling for communication error with picotech

7.18 0.10.1 (2022-01-24)

- **several improvements and fixes for device cube:**
 - privatize `Alarms` and `AlarmsOverview`
 - fix list of cube alarms
 - improve docs
 - fix bugs with earthing sticks
 - fix bug in config dataclass of cube
- introduction of `BoolEnum`
- introduction of `RangeEnum`

- bumpversion -> bump2version

7.19 0.10.0 (2022-01-17)

- Reimplementation of the Cube (before known as Supercube)
- **new names:**
 - Supercube Typ B -> BaseCube
 - Supercube Typ A -> PICube (power inverter Cube)
- **new import:**
 - `from hvl_ccb.dev.supercube import SupercubeB` -> `from hvl_ccb.dev.cube import BaseCube`
- **new programming style:**
 - getter / setter methods -> properties
 - e.g. get: `cube.get_support_output(port=1, contact=1)` -> `cube.support_1.output_1`
 - e.g. set: `cube.get_support_output(port=1, contact=1, state=True)` -> `cube.support_1.output_1 = True`
- unify Exceptions of Cube
- implement Fast Switch-Off of Cube
- remove method `support_output_impulse`
- all active alarms can now be queried `cube.active_alarms()`
- alarms will now result in different logging levels depending on the seriousness of the alarm.
- introduction of limits for slope and safety limit for RedReady
- during the startup the CCB will update the time of the cube.
- verification of inputs
- polarity of DC voltage
- Switch from `python-opcua` to `opcua-asyncio` (former package is no longer maintained)

7.20 0.9.0 (2022-01-07)

- New device: Highland T560 digital delay and pulse generator over Telnet.
- **Rework of the Technix Capacitor Charger.**
 - Moved into a separate sub-package
 - NEW import over `import hvl_ccb.dev.technix as XXX`
 - Slightly adapted behaviour
- Add `validate_tcp_port` to validate port number.
- **Add `validate_and_resolve_host` to validate and resolve host names and IPs.**
 - Remove requirement IPy

- Add a unified CCB Exception schema for all devices and communication protocols.
- Add data conversion functions to README.
- Update CI and devel images from Debian 10 buster to Debian 11 bullseye.
- Fix typing due to numpy update.
- Fix incorrect overloading of `clean_values()` in classes of type `XCommunicationConfig`.

7.21 0.8.5 (2021-11-05)

- Added arbitrary waveform for TiePie signal generation, configurable via `dev.tiepie.generator.TiePieGeneratorConfig.waveform` property.
- In `utils.conversion_sensor`: improvements for class constants; removed SciPy dependency.
- Added Python 3.10 support.

7.22 0.8.4 (2021-10-22)

- `utils.validation.validate_number` extension to handle NumPy arrays and array-like objects.
- `utils.conversion_unit` utility classes handle correctly `NamedTuple` instances.
- `utils.conversion_sensor` and `utils.conversion_unit` code simplification (no `transfer_function_order` attribute) and cleanups.
- Fixed incorrect error logging in `configuration.configdataclass`.
- `comm.telnet.TelnetCommunication` tests fixes for local run errors.

7.23 0.8.3 (2021-09-27)

- New data conversion functions in `utils.conversion_sensor` and `utils.conversion_unit` modules. Note: to use these functions you must install `hvl_ccb` with extra requirement, either `hvl_ccb[conversion]` or `hvl_ccb[all]`.
- Improved documentation with respect to installation of external libraries.

7.24 0.8.2 (2021-08-27)

- **New functionality in `dev.labjack.LabJack`:**
 - configure clock and send timed pulse sequences
 - set DAC/analog output voltage
- Bugfix: ignore random bits sent by to `dev.newport.NewportSMC100PP` controller during start-up/powering-up.

7.25 0.8.1 (2021-08-13)

- Add Python version check (min version error; max version warning).
- Daily checks for upstream dependencies compatibility and devel environment improvements.

7.26 0.8.0 (2021-07-02)

- TCP communication protocol.
- Lauda PRO RP 245 E circulation thermostat device over TCP.
- Pico Technology PT-104 Platinum Resistance Data Logger device as a wrapper of the Python bindings for the PicoSDK.
- In `com.visa.VisaCommunication`: periodic status polling when VISA/TCP keep alive connection is not supported by a host.

7.27 0.7.1 (2021-06-04)

- New `utils.validation` submodule with `validate_bool` and `validate_number` utilities extracted from internal use within a `dev.tiepie` subpackage.
- In `comm.serial.SerialCommunication`:
 - strict encoding errors handling strategy for subclasses,
 - user warning for a low communication timeout value.

7.28 0.7.0 (2021-05-25)

- The `dev.tiepie` module was splitted into a subpackage with, in particular, submodules for each of the device types – `oscilloscope`, `generator`, and `i2c` – and with backward-incompatible direct imports from the submodules.
- In `dev.technix`:
 - fixed communication crash on nested status byte query;
 - added enums for GET and SET register commands.
- Further minor logging improvements: added missing module level logger and removed some error logs in `except` blocks used for a flow control.
- In `examples/` folder renamed consistently all the examples.
- In API documentation: fix incorrect links mapping on inheritance diagrams.

7.29 0.6.1 (2021-05-08)

- In `dev.tiepie`:
 - dynamically set oscilloscope's channel limits in `OscilloscopeChannelParameterLimits`: `input_range` and `trigger_level_abs`, incl. update of latter on each change of `input_range` value of a `TiePieOscilloscopeChannelConfig` instances;
 - quick fix for opening of combined instruments by disabling `OscilloscopeParameterLimits`. `trigger_delay` (an advanced feature);
 - enable automatic devices detection to be able to find network devices with `TiePieOscilloscope`. `list_devices()`.
- Fix `examples/example_labjack.py`.
- Improved logging: consistently use module level loggers, and always log exception tracebacks.
- Improve API documentation: separate pages per modules, each with an inheritance diagram as an overview.

7.30 0.6.0 (2021-04-23)

- Technix capacitor charger using either serial connection or Telnet protocol.
- **Extensions, improvements and fixes in existing devices:**
 - In `dev.tiepie.TiePieOscilloscope`:
 - * redesigned measurement start and data collection API, incl. time out argument, with no/infinite time out option;
 - * trigger allows now a no/infinite time out;
 - * record length and trigger level were fixed to accept, respectively, floating point and integer numbers;
 - * fixed resolution validation bug;
 - `dev.heinzinger.HeinzingerDI` and `dev.rs_rto1024.RTO1024` instances are now resilient to multiple `stop()` calls.
 - In `dev.crylas.CryLasLaser`: default configuration timeout and polling period were adjusted;
 - Fixed PSI9080 example script.
- **Package and source code improvements:**
 - Update to backward-incompatible `pyvisa-py>=0.5.2`. Developers, do update your local development environments!
 - External libraries, like LibTiePie SDK or LJM Library, are now not installed by default; they are now extra installation options.
 - Added Python 3.9 support.
 - Improved number formatting in logs.
 - Typing improvements and fixes for `mypy>=0.800`.

7.31 0.5.0 (2020-11-11)

- TiePie USB oscilloscope, generator and I2C host devices, as a wrapper of the Python bindings for the LibTiePie SDK.
- a FuG Elektronik Power Supply (e.g. Capacitor Charger HCK) using the built-in ADDAT controller with the Probus V protocol over a serial connection
- All devices polling status or measurements use now a `dev.utils.Poller` utility class.
- **Extensions and improvements in existing devices:**
 - In `dev.rs_rto1024.RT01024`: added Channel state, scale, range, position and offset accessors, and measurements activation and read methods.
 - In `dev.sst_luminox.Luminox`: added querying for all measurements in polling mode, and made output mode activation more robust.
 - In `dev.newport.NewportSMC100PP`: an error-prone `wait_until_move_finished` method of replaced by a fixed waiting time, device operations are now robust to a power supply cut, and device restart is not required to apply a start configuration.
- **Other minor improvements:**
 - Single failure-safe starting and stopping of devices sequenced via `dev.base.DeviceSequenceMixin`.
 - Moved `read_text_nonempty` up to `comm.serial.SerialCommunication`.
 - Added development Dockerfile.
 - Updated package and development dependencies: `pymodbus`, `pytest-mock`.

7.32 0.4.0 (2020-07-16)

- **Significantly improved new Supercube device controller:**
 - more robust error-handling,
 - status polling with generic Poller helper,
 - messages and status boards.
 - tested with a physical device,
- Improved OPC UA client wrapper, with better error handling, incl. re-tries on `concurrent.futures.TimeoutError`.
- SST Luminox Oxygen sensor device controller.
- **Backward-incompatible changes:**
 - `CommunicationProtocol.access_lock` has changed type from `threading.Lock` to `threading.RLock`.
 - `ILS2T.relative_step` and `ILS2T.absolute_position` are now called, respectively, `ILS2T.write_relative_step` and `ILS2T.write_absolute_position`.
- **Minor bugfixes and improvements:**
 - fix use of max resolution in `Labjack.set_ain_resolution()`,
 - resolve ILS2T devices relative and absolute position setters race condition,

- added acoustic horn function in the 2015 Supercube.

- **Toolchain changes:**

- add Python 3.8 support,
- drop pytest-runner support,
- ensure compatibility with `labjack_ljm` 2019 version library.

7.33 0.3.5 (2020-02-18)

- Fix issue with reading integers from LabJack LJM Library (device's product ID, serial number etc.)
- Fix development requirements specification (tox version).

7.34 0.3.4 (2019-12-20)

- **New devices using serial connection:**

- Heinzinger Digital Interface I/II and a Heinzinger PNC power supply
- Q-switched Pulsed Laser and a laser attenuator from CryLas
- Newport SMC100PP single axis motion controller for 2-phase stepper motors
- Pfeiffer TPG controller (TPG 25x, TPG 26x and TPG 36x) for Compact pressure Gauges

- PEP 561 compatibility and related corrections for static type checking (now in CI)

- **Refactorings:**

- Protected non-thread safe read and write in communication protocols
- Device sequence mixin: start/stop, add/rm and lookup
- `.format()` to f-strings
- more enumerations and a quite some improvements of existing code

- Improved error docstrings (`:raises:` annotations) and extended tests for errors.

7.35 0.3.3 (2019-05-08)

- Use PyPI labjack-ljm (no external dependencies)

7.36 0.3.2 (2019-05-08)

- INSTALLATION.rst with LJMPython prerequisite info

7.37 0.3.1 (2019-05-02)

- readthedocs.org support

7.38 0.3 (2019-05-02)

- Prevent an automatic close of VISA connection when not used.
- Rhode & Schwarz RTO 1024 oscilloscope using VISA interface over `TCP::INSTR`.
- Extended tests incl. messages sent to devices.
- Added Supercube device using an OPC UA client
- Added Supercube 2015 device using an OPC UA client (for interfacing with old system version)

7.39 0.2.1 (2019-04-01)

- Fix issue with LJMPython not being installed automatically with setuptools.

7.40 0.2.0 (2019-03-31)

- LabJack LJM Library communication wrapper and LabJack device.
- Modbus TCP communication protocol.
- Schneider Electric ILS2T stepper motor drive device.
- Elektro-Automatik PSI9000 current source device and VISA communication wrapper.
- Separate configuration classes for communication protocols and devices.
- Simple experiment manager class.

7.41 0.1.0 (2019-02-06)

- Communication protocol base and serial communication implementation.
- Device base and MBW973 implementation.

**CHAPTER
EIGHT**

INDICES AND TABLES

- genindex
- modindex
- search

PYTHON MODULE INDEX

h

hvl_ccb, 204
hvl_ccb.comm, 35
hvl_ccb.comm.base, 11
hvl_ccb.comm.labjack_ljm, 16
hvl_ccb.comm.modbus_tcp, 19
hvl_ccb.comm.opc, 21
hvl_ccb.comm.serial, 24
hvl_ccb.comm.tcp, 28
hvl_ccb.comm.telnet, 30
hvl_ccb.comm.visa, 32
hvl_ccb.configuration, 200
hvl_ccb.dev, 192
hvl_ccb.dev.base, 187
hvl_ccb.dev.crylas, 46
hvl_ccb.dev.crylas.crylas, 35
hvl_ccb.dev.cube, 61
hvl_ccb.dev.cube.alarms, 46
hvl_ccb.dev.cube.base, 47
hvl_ccb.dev.cube.constants, 53
hvl_ccb.dev.cube.earthing_stick, 55
hvl_ccb.dev.cube.errors, 56
hvl_ccb.dev.cube.picube, 57
hvl_ccb.dev.cube.support, 60
hvl_ccb.dev.cube.switches, 60
hvl_ccb.dev.ea_psi9000, 67
hvl_ccb.dev.ea_psi9000.ea_psi9000, 62
hvl_ccb.dev.fluke884x, 76
hvl_ccb.dev.fluke884x.base, 67
hvl_ccb.dev.fluke884x.constants, 72
hvl_ccb.dev.fluke884x.ranges, 73
hvl_ccb.dev.fug, 89
hvl_ccb.dev.fug.fug, 76
hvl_ccb.dev.heinzinger, 97
hvl_ccb.dev.heinzinger.base, 90
hvl_ccb.dev.heinzinger.constants, 92
hvl_ccb.dev.heinzinger.device, 93
hvl_ccb.dev.heinzinger.mixin, 96
hvl_ccb.dev.highland_t560, 102
hvl_ccb.dev.highland_t560.base, 97
hvl_ccb.dev.highland_t560.channel, 100
hvl_ccb.dev.highland_t560.device, 100

hvl_ccb.dev.labjack, 110
hvl_ccb.dev.labjack.labjack, 103
hvl_ccb.dev.lauda, 118
hvl_ccb.dev.lauda.lauda, 111
hvl_ccb.dev.mbw973, 122
hvl_ccb.dev.mbw973.mbw973, 118
hvl_ccb.dev.newport, 137
hvl_ccb.dev.newport.newport, 123
hvl_ccb.dev.pfeiffer_tpg, 144
hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg, 138
hvl_ccb.dev.protocols, 146
hvl_ccb.dev.protocols.sources, 145
hvl_ccb.dev.rs_rto1024, 154
hvl_ccb.dev.rs_rto1024.rs_rto1024, 146
hvl_ccb.dev.se_ils2t, 162
hvl_ccb.dev.se_ils2t.se_ils2t, 155
hvl_ccb.dev.sst_luminox, 167
hvl_ccb.dev.sst_luminox.sst_luminox, 163
hvl_ccb.dev.technix, 174
hvl_ccb.dev.technix.base, 168
hvl_ccb.dev.technix.device, 171
hvl_ccb.dev.tiepie, 185
hvl_ccb.dev.tiepie.base, 174
hvl_ccb.dev.tiepie.channel, 176
hvl_ccb.dev.tiepie.device, 179
hvl_ccb.dev.tiepie.generator, 180
hvl_ccb.dev.tiepie.oscilloscope, 182
hvl_ccb.dev.tiepie.utils, 185
hvl_ccb.dev.utils, 190
hvl_ccb.dev.visa, 190
hvl_ccb.error, 202
hvl_ccb.experiment_manager, 202
hvl_ccb.utils, 200
hvl_ccb.utils.conversion, 196
hvl_ccb.utils.conversion.map_range, 192
hvl_ccb.utils.conversion.sensor, 193
hvl_ccb.utils.conversion.unit, 194
hvl_ccb.utils.conversion.utils, 195
hvl_ccb.utils.enum, 196
hvl_ccb.utils.poller, 198
hvl_ccb.utils.typing, 199
hvl_ccb.utils.validation, 199

INDEX

A

ABSOLUTE (*TiePieOscilloscopeTriggerLevelMode attribute*), 178
ABSOLUTE_POSITION (*ILS2T.ActionsPtp attribute*), 155
AC (*NewportConfigCommands attribute*), 123
AC_100KV (*PowerSetup attribute*), 54
AC_150KV (*PowerSetup attribute*), 54
AC_200KV (*PowerSetup attribute*), 54
AC_50KV (*PowerSetup attribute*), 54
ac_current_range (*Fluke8845a attribute*), 67
ac_voltage_range (*Fluke8845a attribute*), 67
ACA (*TiePieOscilloscopeChannelCoupling attribute*), 177
acceleration (*NewportSMC100PPConfig attribute*), 131
ACCESS_ENABLE (*ILS2TRegAddr attribute*), 161
access_lock (*CommunicationProtocol attribute*), 14
ACCurrentRange (class in *hvl_ccb.dev.fluke884x.ranges*), 73
ACTION_JOG_VALUE (*ILS2T attribute*), 155
activate_clock_output() (*T560 method*), 100
activate_measurements() (*RTO1024 method*), 146
activate_output() (*Luminox method*), 163
activate_remote_mode() (*Fluke8845a method*), 67
ACTIVE (*CryLasLaser.AnswersStatus attribute*), 39
active_alarms() (*BaseCube method*), 47
ACTIVE_HIGH (*Polarity attribute*), 98
ACTIVE_LOW (*Polarity attribute*), 98
actualsetvalue (*FuGProbusVSetRegisters property*), 85
ACV (*TiePieOscilloscopeChannelCoupling attribute*), 178
ACVoltageRange (class in *hvl_ccb.dev.fluke884x.ranges*), 73
adc_mode (*FuGProbusVMonitorRegisters property*), 85
add_device() (*DeviceSequenceMixin method*), 188
add_device() (*ExperimentManager method*), 202
ADDR_INCORRECT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 133
address (*NewportSMC100PPConfig attribute*), 131
address (*VisaCommunicationConfig property*), 34
address() (*VisaCommunicationConfig.InterfaceType method*), 33

ADMODE (*FuGProbusIVCommands attribute*), 82
ALL (*TiePieOscilloscopeAutoResolutionModes attribute*), 184
all_measurements (*LuminoxMeasurementType attribute*), 165
all_measurements_types() (*LuminoxMeasurementType class method*), 165
amplitude (*TiePieGeneratorConfig property*), 180
ANALOG (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 115
analog_control (*FuGProbusVDIRegisters property*), 84
ANY (*LabJack.DeviceType attribute*), 105
ANY (*LJMCommunicationConfig.ConnectionType attribute*), 17
ANY (*LJMCommunicationConfig.DeviceType attribute*), 17
ANY (*TiePieOscilloscopeTriggerKind attribute*), 178
ApertureRange (class in *hvl_ccb.dev.fluke884x.ranges*), 74
ARBITRARY (*TiePieGeneratorSignalType attribute*), 181
AsyncCommunicationProtocol (class in *hvl_ccb.comm.base*), 11
AsyncCommunicationProtocolConfig (class in *hvl_ccb.comm.base*), 13
ATM (*Pressure attribute*), 194
ATMOSPHERE (*Pressure attribute*), 194
attenuation (*CryLasAttenuator property*), 36
AUTO (*LaudaProRp245eConfig.OperationModeEnum attribute*), 115
AUTO (*RTO1024.TriggerModes attribute*), 146
AUTO (*SwitchOperatingStatus attribute*), 60
auto_install_mode (*T560 property*), 100
auto_install_mode (*T560Config attribute*), 101
auto_laser_on (*CryLasLaserConfig attribute*), 42
auto_resolution_mode (*TiePieOscilloscopeConfig property*), 184
AutoInstallMode (class in *hvl_ccb.dev.highland_t560.base*), 97
AutoNumberNameEnum (class in *hvl_ccb.utils.enum*), 196

B

BA (*NewportConfigCommands* attribute), 123
backlash_compensation (*NewportSMC100PPConfig attribute*), 132
backup_waveform() (*RTO1024* method), 147
BAR (*Pressure* attribute), 194
barometric_pressure (*LuminoxMeasurementType* attribute), 165
base_velocity (*NewportSMC100PPConfig* attribute), 132
BaseCube (*class* in *hvl_ccb.dev(cube.base)*), 47
BaseCubeConfiguration (*class* in *hvl_ccb.dev(cube.base)*), 50
BaseCubeOpcUaCommunication (*class* in *hvl_ccb.dev(cube.base)*), 51
BaseCubeOpcUaCommunicationConfig (*class* in *hvl_ccb.dev(cube.base)*), 51
BATH_TEMP (*LaudaProRp245eCommand* attribute), 113
baudrate (*CryLasAttenuatorSerialCommunicationConfig* attribute), 38
baudrate (*CryLasLaserSerialCommunicationConfig* attribute), 45
baudrate (*FuGSerialCommunicationConfig* attribute), 87
baudrate (*HeinzingerSerialCommunicationConfig* attribute), 91
baudrate (*LuminoxSerialCommunicationConfig* attribute), 166
baudrate (*MBW973SerialCommunicationConfig* attribute), 121
baudrate (*NewportSMC100PPSerialCommunicationConfig* attribute), 136
baudrate (*PfeifferTPGSerialCommunicationConfig* attribute), 143
baudrate (*SerialCommunicationConfig* attribute), 26
BH (*NewportConfigCommands* attribute), 123
board (*VisaCommunicationConfig* attribute), 34
BoolEnum (*class* in *hvl_ccb.utils.enum*), 196
breakdown_detection_active (*BaseCube* property), 47
breakdown_detection_reset() (*BaseCube* method), 47
breakdown_detection_triggered (*BaseCube* property), 47
bufsize (*TcpCommunicationConfig* attribute), 29
build_str() (*LaudaProRp245eCommand* method), 115
BUS (*TriggerSource* attribute), 73
bytesize (*CryLasAttenuatorSerialCommunicationConfig* attribute), 38
bytesize (*CryLasLaserSerialCommunicationConfig* attribute), 45
bytesize (*FuGSerialCommunicationConfig* attribute), 87

bytesize (*HeinzingerSerialCommunicationConfig* attribute), 91
bytesize (*LuminoxSerialCommunicationConfig* attribute), 166
bytesize (*MBW973SerialCommunicationConfig* attribute), 121
bytesize (*NewportSMC100PPSerialCommunicationConfig* attribute), 136
bytesize (*PfeifferTPGSerialCommunicationConfig* attribute), 143
Bytesize (*SerialCommunicationConfig* attribute), 26
bytesize (*SerialCommunicationConfig* attribute), 26
C
C (*LabJack.TemperatureUnit* attribute), 105
C (*LabJack.ThermocoupleType* attribute), 106
C (*Temperature* attribute), 195
calibration_factor (*CryLasLaserConfig* attribute), 42
calibration_factor (*LEM4000S* attribute), 193
calibration_mode (*FuGProbusVDIRegisters* property), 84
cc_mode (*FuGProbusVDIRegisters* property), 84
CCBError, 202
cee16_socket (*BaseCube* property), 47
CELSIUS (*Temperature* attribute), 195
ch_a (*T560* property), 100
ch_b (*T560* property), 100
ch_c (*T560* property), 101
ch_d (*T560* property), 101
channels_enabled (*TiePieOscilloscope* property), 182
check_for_error() (*NewportSMC100PPSerialCommunication* method), 134
check_master_slave_config() (*PSI9000* method), 62
chunk_size (*VisaCommunicationConfig* attribute), 34
clean_amplitude() (*TiePieGeneratorConfig* method), 180
clean_auto_resolution_mode() (*TiePieOscilloscopeConfig* static method), 184
clean_coupling() (*TiePieOscilloscopeChannelConfig* static method), 177
clean_enabled() (*TiePieGeneratorConfig* static method), 180
clean_enabled() (*TiePieOscilloscopeChannelConfig* static method), 177
clean_frequency() (*TiePieGeneratorConfig* method), 180
clean_input_range() (*TiePieOscilloscopeChannelConfig* method), 177
clean_offset() (*TiePieGeneratorConfig* method), 180
clean_pre_sample_ratio() (*TiePieOscilloscopeConfig* method), 184

clean_probe_offset() (*TiePieOscilloscopeChannelConfig method*), 177
clean_record_length() (*TiePieOscilloscopeConfig method*), 184
clean_resolution() (*TiePieOscilloscopeConfig static method*), 184
clean_sample_rate() (*TiePieOscilloscopeConfig method*), 184
clean_signal_type() (*TiePieGeneratorConfig static method*), 180
clean_trigger_enabled() (*TiePieOscilloscopeChannelConfig static method*), 177
clean_trigger_hysteresis() (*TiePieOscilloscopeChannelConfig method*), 177
clean_trigger_kind() (*TiePieOscilloscopeChannelConfig static method*), 177
clean_trigger_level() (*TiePieOscilloscopeChannelConfig method*), 177
clean_trigger_level_mode() (*TiePieOscilloscopeChannelConfig static method*), 177
clean_trigger_timeout() (*TiePieOscilloscopeConfig method*), 184
clean_values() (*AsyncCommunicationProtocolConfig method*), 13
clean_values() (*BaseCubeConfiguration method*), 50
clean_values() (*CryLasAttenuatorConfig method*), 36
clean_values() (*CryLasLaserConfig method*), 42
clean_values() (*EmptyConfig method*), 189, 200
clean_values() (*Fluke8845aConfig method*), 69
clean_values() (*FuGConfig method*), 78
clean_values() (*HeinzingerConfig method*), 95
clean_values() (*ILS2TConfig method*), 159
clean_values() (*LaudaProRp245eConfig method*), 115
clean_values() (*LaudaProRp245eTcpCommunicationConfig method*), 117
clean_values() (*LJMCommunicationConfig method*), 18
clean_values() (*LuminoxConfig method*), 164
clean_values() (*MBW973Config method*), 120
clean_values() (*ModbusTcpCommunicationConfig method*), 20
clean_values() (*NewportSMC100PPConfig method*), 132
clean_values() (*OpcUaCommunicationConfig method*), 22
clean_values() (*PfeifferTPGConfig method*), 141
clean_values() (*PICubeConfiguration method*), 58
clean_values() (*PSI9000Config method*), 64
clean_values() (*SerialCommunicationConfig method*), 26
clean_values() (*T560Config method*), 101
clean_values() (*TcpCommunicationConfig method*), 29
clean_values() (*TechnixConfig method*), 172
clean_values() (*TelnetCommunicationConfig method*), 31
clean_values() (*TiePieDeviceConfig method*), 174
clean_values() (*VisaCommunicationConfig method*), 34
clean_waveform() (*TiePieGeneratorConfig method*), 180
clear_display_message() (*Fluke8845a method*), 67
clear_error_queue() (*Fluke8845a method*), 68
Client (class in *hvl_ccb.comm.opc*), 21
CLOSE (*SwitchOperation attribute*), 61
close() (*CommunicationProtocol method*), 14
close() (*LaudaProRp245eTcpCommunication method*), 117
close() (*LJMCommunication method*), 16
close() (*ModbusTcpCommunication method*), 19
close() (*NullCommunicationProtocol method*), 14
close() (*OpcUaCommunication method*), 21
close() (*SerialCommunication method*), 24
close() (*Tcp method*), 28
close() (*TelnetCommunication method*), 30
close() (*VisaCommunication method*), 32
close_shutter() (*CryLasLaser method*), 40
CLOSED (*CryLasLaser.AnswersShutter attribute*), 39
CLOSED (*CryLasLaserShutterStatus attribute*), 46
CLOSED (*DoorStatus attribute*), 53
CLOSED (*SwitchStatus attribute*), 61
CMD_EXEC_ERROR (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_CC (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_CONFIGURATION (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_DISABLE (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_HOMING (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_MOVING (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_NOT_REFERENCED (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
CMD_NOT_ALLOWED_PP (*New-*

`portSMC100PPSerialCommunication.ControllerErrors` method), 160
attribute), 133
`CMD_NOT_ALLOWED_READY` (New-
`portSMC100PPSerialCommunication.ControllerErrors` static method), 117
attribute), 133
`CMR (PfeifferTPG.SensorTypes attribute)`, 139
`CODE_OR_ADDR_INVALID` (New-
`portSMC100PPSerialCommunication.ControllerErrors` method), 166
attribute), 133
`collect_measurement_data()` (`TiePieOscilloscope`
method), 182
`com (SingleCommDevice property)`, 189
`COM_TIME_OUT` (`LaudaProRp245eCommand` attribute),
114
`com_time_out` (`LaudaProRp245eConfig` attribute), 115
`COM_TIMEOUT` (`NewportSMC100PPSerialCommunication.ControllerErrors` method), 134
attribute), 133
`command (LuminoxMeasurementType property)`, 165
`COMMAND (TriggerMode attribute)`, 99
`command()` (`FuGProbusIV` method), 81
`communication_channel` (`TechnixConfig` attribute),
172
`CommunicationError`, 14
`CommunicationProtocol` (class
`hvl_ccb.comm.base`), 14
`config (ConfigurationMixin property)`, 200
`CONFIG (FuGProbusVRegisterGroups attribute)`, 85
`CONFIG (NewportSMC100PP.StateMessages attribute)`,
124
`CONFIG (NewportStates attribute)`, 137
`config_cls()` (`AsyncCommunicationProtocol` static
method), 11
`config_cls()` (`BaseCube` static method), 48
`config_cls()` (`BaseCubeOpcUaCommunication` static
method), 51
`config_cls()` (`ConfigurationMixin` static method), 200
`config_cls()` (`CryLasAttenuator` static method), 36
`config_cls()` (`CryLasAttenuatorSerialCommunication`
static method), 37
`config_cls()` (`CryLasLaser` static method), 40
`config_cls()` (`CryLasLaserSerialCommunication`
static method), 44
`config_cls()` (`Device` static method), 187
`config_cls()` (`Fluke8845a` static method), 68
`config_cls()` (`Fluke8845aTelnetCommunication` static
method), 70
`config_cls()` (`FuGProbusIV` static method), 81
`config_cls()` (`FuGSerialCommunication` static
method), 87
`config_cls()` (`Heinzinger` static method), 94
`config_cls()` (`HeinzingerSerialCommunication` static
method), 90
`config_cls()` (`ILS2T` static method), 156
`config_cls()` (`ILS2TModbusTcpCommunication` static
method), 160
`config_cls()` (`LaudaProRp245e` static method), 111
`config_cls()` (`LaudaProRp245eTcpCommunication`
static method), 117
`config_cls()` (`LJMCommunication` static method), 16
`config_cls()` (`Luminox` static method), 163
`config_cls()` (`LuminoxSerialCommunication` static
method), 121
`config_cls()` (`ModbusTcpCommunication` static
method), 19
`config_cls()` (`NewportSMC100PP` static method), 125
`config_cls()` (`NewportSMC100PPSerialCommunication`
static method), 134
`config_cls()` (`NullCommunicationProtocol` static
method), 14
`config_cls()` (`OpcUaCommunication` static method),
21
`config_cls()` (`PfeifferTPG` static method), 139
`config_cls()` (`PfeifferTPGSerialCommunication` static
method), 142
`config_cls()` (`PICube` static method), 57
`config_cls()` (`PICubeOpcUaCommunication` static
method), 59
`config_cls()` (`PSI9000` static method), 62
`config_cls()` (`PSI9000VisaCommunication` static
method), 66
`config_cls()` (`RTO1024` static method), 147
`config_cls()` (`RTO1024VisaCommunication` static
method), 153
`config_cls()` (`SerialCommunication` static method), 25
`config_cls()` (`SyncCommunicationProtocol` static
method), 15
`config_cls()` (`T560` static method), 101
`config_cls()` (`T560Communication` static method), 98
`config_cls()` (`Tcp` static method), 28
`config_cls()` (`Technix` static method), 171
`config_cls()` (`TechnixSerialCommunication` static
method), 168
`config_cls()` (`TechnixTelnetCommunication` static
method), 169
`config_cls()` (`TelnetCommunication` static method), 30
`config_cls()` (`TiePieOscilloscope` static method), 182
`config_cls()` (`VisaCommunication` static method), 32
`config_cls()` (`VisaDevice` static method), 190
`config_gen` (`TiePieGeneratorMixin` attribute), 181
`config_high_pulse()` (`LabJack` method), 106
`config_osc` (`TiePieOscilloscope` attribute), 182
`config_osc_channel_dict` (`TiePieOscilloscope`
attribute), 183
`config_status` (`FuG` property), 76
`configdataclass()` (in module `hvl_ccb.configuration`),

201
configuration_save_json() (*ConfigurationMixin method*), 200
ConfigurationMixin (*class in hvl_ccb.configuration*), 200
ConfigurationValueWarning, 200
connection_type (*LJMCommunicationConfig attribute*), 18
CONT_MODE (*LaudaProRp245eCommand attribute*), 114
continue_ramp() (*LaudaProRp245e method*), 111
control_mode (*LaudaProRp245eConfig attribute*), 116
CONVERSION (*LEM4000S attribute*), 193
convert() (*LEM4000S method*), 193
convert() (*LMT70A method*), 194
convert() (*Sensor method*), 194
convert() (*Unit class method*), 195
convert_to_range1() (*MapRanges method*), 193
convert_to_range2() (*MapRanges method*), 193
convert_value_to_str() (*in module hvl_ccb.utils.conversion.utils*), 195
ConvertibleTypes (*in module hvl_ccb.utils.typing*), 199
COOLOFF (*LaudaProRp245eConfig.OperationModeEnum attribute*), 115
COOLON (*LaudaProRp245eConfig.OperationModeEnum attribute*), 115
coupling (*TiePieOscilloscopeChannelConfig property*), 177
CR (*FuGTerminators attribute*), 88
create_serial_port() (*SerialCommunicationConfig method*), 26
create_telnet() (*TelnetCommunicationConfig method*), 31
CRLF (*FuGTerminators attribute*), 88
CryLasAttenuator (*class in hvl_ccb.dev.crylas.crylas*), 36
CryLasAttenuatorConfig (*class in hvl_ccb.dev.crylas.crylas*), 36
CryLasAttenuatorError, 37
CryLasAttenuatorSerialCommunication (*class in hvl_ccb.dev.crylas.crylas*), 37
CryLasAttenuatorSerialCommunicationConfig (*class in hvl_ccb.dev.crylas.crylas*), 37
CryLasLaser (*class in hvl_ccb.dev.crylas.crylas*), 39
CryLasLaser.AnswersShutter (*class in hvl_ccb.dev.crylas.crylas*), 39
CryLasLaser.AnswersStatus (*class in hvl_ccb.dev.crylas.crylas*), 39
CryLasLaser.LaserStatus (*class in hvl_ccb.dev.crylas.crylas*), 39
CryLasLaser.RepetitionRates (*class in hvl_ccb.dev.crylas.crylas*), 39
CryLasLaserConfig (*class in hvl_ccb.dev.crylas.crylas*), 42
CryLasLaserError, 43
CryLasLaserNotReadyError, 43
CryLasLaserPoller (*class in hvl_ccb.dev.crylas.crylas*), 43
CryLasLaserSerialCommunication (*class in hvl_ccb.dev.crylas.crylas*), 43
CryLasLaserSerialCommunicationConfig (*class in hvl_ccb.dev.crylas.crylas*), 44
CryLasLaserShutterStatus (*class in hvl_ccb.dev.crylas.crylas*), 46
CubeEarthingStickOperationError, 56
CubeError, 56
CubeRemoteControlError, 56
CubeStatusChangeError, 56
CubeStopError, 56
current (*FuG property*), 76
CURRENT (*FuGProbusIVCommands attribute*), 82
CURRENT (*FuGReadbackChannels attribute*), 87
current (*Heinzinger property*), 94
current (*Source property*), 145
current (*Technix property*), 171
CURRENT_AC (*MeasurementFunction attribute*), 72
CURRENT_DC (*MeasurementFunction attribute*), 72
current_filter (*Fluke8845a attribute*), 68
current_lower_limit (*PSI9000Config attribute*), 65
current_monitor (*FuG property*), 76
current_primary (*PICube property*), 57
current_upper_limit (*PSI9000Config attribute*), 65
CurrentPosition (*New-portSMC100PPConfig.HomeSearch attribute*), 131
cv_mode (*FuGProbusVDIRegisters property*), 84

D

datachange_notification() (*OpcUaSubHandler method*), 24
date_of_manufacture (*LuminoxMeasurementType attribute*), 165
datetime_to_opc() (*BaseCube class method*), 48
DC (*TiePieGeneratorSignalType attribute*), 181
DC_140KV (*PowerSetup attribute*), 55
DC_280KV (*PowerSetup attribute*), 55
dc_current_range (*Fluke8845a attribute*), 68
dc_voltage_range (*Fluke8845a attribute*), 68
DC_VOLTAGE_TOO_LOW (*New-portSMC100PP.MotorErrors attribute*), 124
DCA (*TiePieOscilloscopeChannelCoupling attribute*), 178
DCCurrentRange (*class in hvl_ccb.dev.fluke884x.ranges*), 74
DCV (*TiePieOscilloscopeChannelCoupling attribute*), 178
DCVoltageRange (*class in hvl_ccb.dev.fluke884x.ranges*), 74
default_com_cls() (*BaseCube static method*), 48

default_com_cls() (*CryLasAttenuator static method*), 36
default_com_cls() (*CryLasLaser static method*), 40
default_com_cls() (*Fluke8845a static method*), 68
default_com_cls() (*FuGProbusIV static method*), 81
default_com_cls() (*Heinzinger static method*), 94
default_com_cls() (*ILS2T static method*), 156
default_com_cls() (*LabJack static method*), 106
default_com_cls() (*LaudaProRp245e static method*), 111
default_com_cls() (*Luminox static method*), 163
default_com_cls() (*MBW973 static method*), 119
default_com_cls() (*NewportSMC100PP static method*), 125
default_com_cls() (*PfeifferTPG static method*), 139
default_com_cls() (*PICube static method*), 57
default_com_cls() (*PSI9000 static method*), 62
default_com_cls() (*RTO1024 static method*), 147
default_com_cls() (*SingleCommDevice static method*), 189
default_com_cls() (*T560 static method*), 101
default_com_cls() (*Technix method*), 171
default_com_cls() (*TiePieOscilloscope static method*), 183
default_com_cls() (*VisaDevice static method*), 190
DEFAULT_IO_SCANNING_CONTROL_VALUES (*ILS2T attribute*), 155
default_n_attempts_read_text_nonempty (*AsyncCommunicationProtocolConfig attribute*), 13
default_n_attempts_read_text_nonempty (*FuGSerialCommunicationConfig attribute*), 88
default_n_attempts_read_text_nonempty (*HeinzingerSerialCommunicationConfig attribute*), 91
default_number_of_recordings (*HeinzingerConfig attribute*), 95
DeprecatedHeinzingerMixin (class in *hvl_ccb.dev.heinzinger.mixin*), 96
Device (class in *hvl_ccb.dev.base*), 187
DEVICE_TYPE (*LaudaProRp245eCommand attribute*), 114
device_type (*LJMCommunicationConfig attribute*), 18
DeviceError, 187
DeviceExistingError, 187
DeviceFailuresError, 187
devices_failed_start (*DeviceSequenceMixin attribute*), 188
devices_failed_start (*ExperimentManager attribute*), 203
devices_failed_stop (*DeviceSequenceMixin attribute*), 188
devices_failed_stop (*ExperimentManager attribute*), 203
DeviceSequenceMixin (class in *hvl_ccb.dev.base*), 187
di (*FuG property*), 77
digital_control (*FuGProbusVDIRegisters property*), 84
DIOChannel (*LabJack attribute*), 105
DIODE (*MeasurementFunction attribute*), 72
DISABLE (*NewportStates attribute*), 137
disable() (*ILS2T method*), 156
DISABLE_FROM_JOGGING (*portSMC100PP.StateMessages attribute*), 124
DISABLE_FROM_MOVING (*portSMC100PP.StateMessages attribute*), 124
DISABLE_FROM_READY (*portSMC100PP.StateMessages attribute*), 125
disable_pulses() (*LabJack method*), 106
DISABLED (*TiePieOscilloscopeAutoResolutionModes attribute*), 184
DisableEspStageCheck (*portSMC100PPConfig.EspStageConfig attribute*), 131
disarm_trigger() (*T560 method*), 101
disconnect() (*Client method*), 21
DISPLACEMENT_OUT_OF_LIMIT (*portSMC100PPSerialCommunication.ControllerErrors attribute*), 133
display_enable (*Fluke8845a property*), 68
DISPLAY_MAX_LENGTH (*Fluke8845a attribute*), 67
display_message (*Fluke8845a property*), 68
display_message_board() (*BaseCube method*), 48
display_status_board() (*BaseCube method*), 48
do_ioscanning_write() (*ILS2T method*), 157
door_1_status (*BaseCube attribute*), 48
door_2_status (*BaseCube attribute*), 48
door_3_status (*BaseCube attribute*), 48
DoorStatus (class in *hvl_ccb.dev.constants*), 53

E

E (*LabJack.ThermocoupleType attribute*), 106
E0 (*FuGErrorcodes attribute*), 79
E1 (*FuGErrorcodes attribute*), 79
E10 (*FuGErrorcodes attribute*), 79
E100 (*FuGErrorcodes attribute*), 79
E106 (*FuGErrorcodes attribute*), 79
E11 (*FuGErrorcodes attribute*), 79
E115 (*FuGErrorcodes attribute*), 79
E12 (*FuGErrorcodes attribute*), 79
E125 (*FuGErrorcodes attribute*), 79
E13 (*FuGErrorcodes attribute*), 79
E135 (*FuGErrorcodes attribute*), 79
E14 (*FuGErrorcodes attribute*), 80
E145 (*FuGErrorcodes attribute*), 80
E15 (*FuGErrorcodes attribute*), 80

E155 (*FuGErrorcodes attribute*), 80
E16 (*FuGErrorcodes attribute*), 80
E165 (*FuGErrorcodes attribute*), 80
E2 (*FuGErrorcodes attribute*), 80
E206 (*FuGErrorcodes attribute*), 80
E306 (*FuGErrorcodes attribute*), 80
E4 (*FuGErrorcodes attribute*), 80
E5 (*FuGErrorcodes attribute*), 80
E504 (*FuGErrorcodes attribute*), 80
E505 (*FuGErrorcodes attribute*), 80
E6 (*FuGErrorcodes attribute*), 80
E666 (*FuGErrorcodes attribute*), 80
E7 (*FuGErrorcodes attribute*), 80
E8 (*FuGErrorcodes attribute*), 80
E9 (*FuGErrorcodes attribute*), 80
earthing_rod_1_status (*BaseCube attribute*), 48
earthing_rod_2_status (*BaseCube attribute*), 48
earthing_rod_3_status (*BaseCube attribute*), 48
EarthingRodStatus (class in *hvl_ccb.dev(cube.constants)*), 53
EEPROM_ACCESS_ERROR (New-
portSMC100PPSerialCommunication.ControllerErrors attribute), 134
EIGHT (*RecordingsEnum attribute*), 93
EIGHT_BIT (*TiePieOscilloscopeResolution attribute*), 185
EIGHT_HUNDRED_MILLI_VOLT (*TiePieOscilloscop eRange attribute*), 178
EIGHT_VOLT (*TiePieOscilloscopeRange attribute*), 178
EIGHTBITS (*SerialCommunicationBytesize attribute*), 25
EIGHTY_VOLT (*TiePieOscilloscopeRange attribute*), 178
EmptyConfig (class in *hvl_ccb.configuration*), 200
EmptyConfig (class in *hvl_ccb.dev.base*), 189
enable() (*ILS2T method*), 157
enable_clock() (*LabJack method*), 106
enabled (*TiePieGeneratorConfig property*), 180
enabled (*TiePieOscilloscopeChannelConfig property*), 177
EnableEspStageCheck (New-
portSMC100PPConfig.EspStageConfig attribute), 131
encoding (*AsyncCommunicationProtocolConfig attribute*), 13
encoding (*NewportSMC100PPSerialCommunicationConfig attribute*), 136
encoding_error_handling (*AsyncCommunicationProtocolConfig attribute*), 13
encoding_error_handling (New-
portSMC100PPSerialCommunicationConfig attribute), 136
EndOfRunSwitch (New-
portSMC100PPConfig.HomeSearch attribute), 131
EndOfRunSwitch_and_Index (New-
portSMC100PPConfig.HomeSearch attribute), 131
portSMC100PPConfig.HomeSearch attribute), 131
endpoint_name (*BaseCubeOpcUaCommunicationConfig attribute*), 52
endpoint_name (*OpcUaCommunicationConfig attribute*), 23
endpoint_name (*PICubeOpcUaCommunicationConfig attribute*), 59
ERROR (*DoorStatus attribute*), 53
ERROR (*ExperimentStatus attribute*), 203
ERROR (*SafetyStatus attribute*), 55
ERROR (*SwitchStatus attribute*), 61
errorcode (*FuGError attribute*), 79
ESP_STAGE_NAME_INVALID (New-
portSMC100PPSerialCommunication.ControllerErrors attribute), 134
ETH (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 115
in ETHERNET (*LJMCommunicationConfig.ConnectionType attribute*), 17
EVEN (*SerialCommunicationParity attribute*), 27
Event_notification() (*OpcUaSubHandler method*), 24
EXECUTE (*FuGProbusIVCommands attribute*), 82
execute_absolute_position() (*ILS2T method*), 157
execute_on_x (*FuGProbusVConfigRegisters property*), 83
execute_relative_step() (*ILS2T method*), 157
EXECUTEONX (*FuGProbusIVCommands attribute*), 82
exit_configuration() (*NewportSMC100PP method*), 125
exit_configuration_wait_sec (New-
portSMC100PPConfig attribute), 132
EXPERIMENT_BLOCKED (*EarthingRodStatus attribute*), 53
EXPERIMENT_READY (*EarthingRodStatus attribute*), 53
ExperimentError, 202
ExperimentManager (class in *hvl_ccb.experiment_manager*), 202
ExperimentStatus (class in *hvl_ccb.experiment_manager*), 203
EXPT100 (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 115
EXT_FALLING_EDGE (*TriggerMode attribute*), 99
EXT_RISING_EDGE (*TriggerMode attribute*), 99
EXTERNAL (*TriggerSource attribute*), 73
EXTERNAL_SOURCE (*PowerSetup attribute*), 55
EXTERNAL_TEMP (*LaudaProRp245eCommand attribute*), 114
F
F (*LabJack.TemperatureUnit attribute*), 105
F (*Temperature attribute*), 195
FAHRENHEIT (*Temperature attribute*), 195
failures (*DeviceFailuresError attribute*), 187

FALLING (*TiePieOscilloscopeTriggerKind* attribute), 178
FAST (*ILS2T.Ref16Jog* attribute), 156
FAST_FILTER (*FilterRange* attribute), 75
fetch() (*Fluke8845a* method), 68
file_copy() (*RTO1024* method), 147
FilterRange (class in *hvl_ccb.dev.fluke884x.ranges*), 75
finish() (*ExperimentManager* method), 203
FINISHED (*ExperimentStatus* attribute), 203
FINISHING (*ExperimentStatus* attribute), 203
fire_trigger() (*T560* method), 101
FIRMWARE (*FuGRreadbackChannels* attribute), 87
FIVE_MHZ (*LabJack.ClockFrequency* attribute), 104
FIVEBITS (*SerialCommunicationBytesize* attribute), 25
FLT_INFO (*ILS2TRegAddr* attribute), 161
FLT_MEM_DEL (*ILS2TRegAddr* attribute), 161
FLT_MEM_RESET (*ILS2TRegAddr* attribute), 161
Fluke8845a (class in *hvl_ccb.dev.fluke884x.base*), 67
Fluke8845aCheckError, 72
Fluke8845aConfig (class in *hvl_ccb.dev.fluke884x.base*), 69
Fluke8845aError, 72
Fluke8845aTelnetCommunication (class in *hvl_ccb.dev.fluke884x.base*), 70
Fluke8845aTelnetCommunicationConfig (class in *hvl_ccb.dev.fluke884x.base*), 70
Fluke8845aUnknownCommandError, 72
FOLLOWING_ERROR (*NewportSMC100PP.MotorErrors* attribute), 124
FOLLOWRAMP (*FuGRampModes* attribute), 86
force_trigger() (*TiePieOscilloscope* method), 183
force_value() (*AsyncCommunicationProtocolConfig* method), 13
force_value() (*BaseCubeConfiguration* method), 50
force_value() (*BaseCubeOpcUaCommunicationConfig* method), 52
force_value() (*CryLasAttenuatorConfig* method), 36
force_value() (*CryLasAttenuatorSerialCommunicationConfig* method), 38
force_value() (*CryLasLaserConfig* method), 42
force_value() (*CryLasLaserSerialCommunicationConfig* method), 45
force_value() (*EmptyConfig* method), 189, 201
force_value() (*Fluke8845aConfig* method), 69
force_value() (*Fluke8845aTelnetCommunicationConfig* method), 71
force_value() (*FuGConfig* method), 78
force_value() (*FuGSerialCommunicationConfig* method), 88
force_value() (*HeinzingerConfig* method), 95
force_value() (*HeinzingerSerialCommunicationConfig* method), 91
force_value() (*ILS2TConfig* method), 159
force_value() (*ILS2TModbusTcpCommunicationConfig* method), 160
force_value() (*LaudaProRp245eConfig* method), 116
force_value() (*LaudaProRp245eTcpCommunicationConfig* method), 117
force_value() (*LJMCommunicationConfig* method), 18
force_value() (*LuminoxConfig* method), 164
force_value() (*LuminoxSerialCommunicationConfig* method), 166
force_value() (*MBW973Config* method), 120
force_value() (*MBW973SerialCommunicationConfig* method), 121
force_value() (*ModbusTcpCommunicationConfig* method), 20
force_value() (*NewportSMC100PPConfig* method), 132
force_value() (*NewportSMC100PPSerialCommunicationConfig* method), 136
force_value() (*OpcUaCommunicationConfig* method), 23
force_value() (*PfeifferTPGConfig* method), 141
force_value() (*PfeifferTPGSerialCommunicationConfig* method), 143
force_value() (*PICubeConfiguration* method), 58
force_value() (*PICubeOpcUaCommunicationConfig* method), 59
force_value() (*PSI9000Config* method), 65
force_value() (*PSI9000VisaCommunicationConfig* method), 66
force_value() (*RTO1024Config* method), 153
force_value() (*RTO1024VisaCommunicationConfig* method), 154
force_value() (*SerialCommunicationConfig* method), 26
force_value() (*T560CommunicationConfig* method), 99
force_value() (*T560Config* method), 102
force_value() (*TcpCommunicationConfig* method), 29
force_value() (*TechnixConfig* method), 172
force_value() (*TechnixSerialCommunicationConfig* method), 169
force_value() (*TechnixTelnetCommunicationConfig* method), 170
force_value() (*TelnetCommunicationConfig* method), 31
force_value() (*TiePieDeviceConfig* method), 175
force_value() (*VisaCommunicationConfig* method), 34
force_value() (*VisaDeviceConfig* method), 191
FORTY_MHZ (*LabJack.ClockFrequency* attribute), 104
FORTY_VOLT (*TiePieOscilloscopeRange* attribute), 178
FOUR (*RecordingsEnum* attribute), 93
FOUR_HUNDRED_MILLI_AMPERE (*ACCurrantRange* attribute), 73

FOUR_HUNDRED_MILLI_AMPERE (<i>DCCurrentRange attribute</i>), 74		gate_polarity (<i>T560 property</i>), 101
FOUR_HUNDRED_MILLI_VOLT (<i>TiePieOscilloscopeRange attribute</i>), 178		GateMode (<i>class in hvl_ccb.dev.highland_t560.base</i>), 98
FOUR_VOLT (<i>TiePieOscilloscopeRange attribute</i>), 178		GENERATOR (<i>TiePieDeviceType attribute</i>), 175
FOUR_WIRE_RESISTANCE (<i>MeasurementFunction attribute</i>), 72		generator_start() (<i>TiePieGeneratorMixin method</i>), 181
four_wire_resistance_range (<i>Fluke8845a attribute</i>), 68		generator_stop() (<i>TiePieGeneratorMixin method</i>), 181
FOURTEEN_BIT (<i>TiePieOscilloscopeResolution attribute</i>), 185		get_acceleration() (<i>NewportSMC100PP method</i>), 126
FREERUN (<i>RTO1024.TriggerModes attribute</i>), 146		get_acquire_length() (<i>RTO1024 method</i>), 147
FREQUENCY (<i>MeasurementFunction attribute</i>), 72		get_ain() (<i>LabJack method</i>), 106
frequency (<i>PICube property</i>), 57		get_bath_temp() (<i>LaudaProRp245e method</i>), 111
frequency (<i>T560 property</i>), 101		get_by_p_id() (<i>LabJack.DeviceType class method</i>), 105
frequency (<i>TiePieGeneratorConfig property</i>), 180		get_by_p_id() (<i>LJMCommunicationConfig.DeviceType class method</i>), 17
frequency_aperture (<i>Fluke8845a attribute</i>), 68		get_cal_current_source() (<i>LabJack method</i>), 107
FRM (<i>NewportConfigCommands attribute</i>), 123		get_channel_offset() (<i>RTO1024 method</i>), 147
from_json() (<i>ConfigurationMixin class method</i>), 200		get_channel_position() (<i>RTO1024 method</i>), 148
FRS (<i>NewportConfigCommands attribute</i>), 123		get_channel_range() (<i>RTO1024 method</i>), 148
FuG (<i>class in hvl_ccb.dev.fug.fug</i>), 76		get_channel_scale() (<i>RTO1024 method</i>), 148
FuGConfig (<i>class in hvl_ccb.dev.fug.fug</i>), 78		get_channel_state() (<i>RTO1024 method</i>), 148
FuGDigitalVal (<i>class in hvl_ccb.dev.fug.fug</i>), 79		get_clock() (<i>LabJack method</i>), 107
FuGError, 79		get_controller_information() (<i>NewportportSMC100PP method</i>), 126
FuGErrorcodes (<i>class in hvl_ccb.dev.fug.fug</i>), 79		get_dc_volt() (<i>ILS2T method</i>), 157
FuGMonitorModes (<i>class in hvl_ccb.dev.fug.fug</i>), 80		get_device() (<i>DeviceSequenceMixin method</i>), 188
FuGPolarities (<i>class in hvl_ccb.dev.fug.fug</i>), 81		get_device_by_serial_number() (<i>in module hvl_ccb.dev.tiepie.base</i>), 175
FuGProbusIV (<i>class in hvl_ccb.dev.fug.fug</i>), 81		get_device_type() (<i>LaudaProRp245e method</i>), 111
FuGProbusIVCommands (<i>class in hvl_ccb.dev.fug.fug</i>), 82		get_devices() (<i>DeviceSequenceMixin method</i>), 188
FuGProbusV (<i>class in hvl_ccb.dev.fug.fug</i>), 82		get_digital_input() (<i>LabJack method</i>), 107
FuGProbusVConfigRegisters (<i>class in hvl_ccb.dev.fug.fug</i>), 83	in	get_error_code() (<i>ILS2T method</i>), 157
FuGProbusVDIRegisters (<i>class in hvl_ccb.dev.fug.fug</i>), 84		get_error_queue() (<i>VisaDevice method</i>), 191
FuGProbusVDORRegisters (<i>class in hvl_ccb.dev.fug.fug</i>), 84		get_full_scale_mbar() (<i>PfeifferTPG method</i>), 139
FuGProbusVMonitorRegisters (<i>class in hvl_ccb.dev.fug.fug</i>), 85	in	get_full_scale_unitless() (<i>PfeifferTPG method</i>), 139
FuGProbusVRegisterGroups (<i>class in hvl_ccb.dev.fug.fug</i>), 85	in	get_identification() (<i>VisaDevice method</i>), 191
FuGProbusVSetRegisters (<i>class in hvl_ccb.dev.fug.fug</i>), 85	in	get_interface_version() (<i>Heinzinger method</i>), 94
FuGRampModes (<i>class in hvl_ccb.dev.fug.fug</i>), 86		get_motor_configuration() (<i>NewportSMC100PP method</i>), 126
FuGReadbackChannels (<i>class in hvl_ccb.dev.fug.fug</i>), 86		get_move_duration() (<i>NewportSMC100PP method</i>), 126
FuGSerialCommunication (<i>class in hvl_ccb.dev.fug.fug</i>), 87	in	get_negative_software_limit() (<i>NewportportSMC100PP method</i>), 127
FuGSerialCommunicationConfig (<i>class in hvl_ccb.dev.fug.fug</i>), 87	in	get_number_of_recordings() (<i>Deprecated-HeinzingerMixin static method</i>), 96
FuGTerminators (<i>class in hvl_ccb.dev.fug.fug</i>), 88		get_objects_node() (<i>Client method</i>), 21
G		get_objects_node() (<i>Server method</i>), 24
gate_mode (<i>T560 property</i>), 101		get_output() (<i>PSI9000 method</i>), 63
		get_position() (<i>ILS2T method</i>), 157
		get_position() (<i>NewportSMC100PP method</i>), 127
		get_positive_software_limit() (<i>NewportportSMC100PP method</i>), 127

get_product_id() (*LabJack method*), 107
 get_product_name() (*LabJack method*), 107
 get_product_type() (*LabJack method*), 107
 get_pulse_energy_and_rate() (*CryLasLaser method*), 40
 get_reference_point() (*RTO1024 method*), 148
 get_register() (*FuGProbusV method*), 82
 get_repetitions() (*RTO1024 method*), 148
 get_sbus_rh() (*LabJack method*), 107
 get_sbus_temp() (*LabJack method*), 108
 get_serial_number() (*Heinzinger method*), 94
 get_serial_number() (*LabJack method*), 108
 get_state() (*NewportSMC100PP method*), 127
 get_status() (*ILS2T method*), 157
 get_system_lock() (*PSI9000 method*), 63
 get_temperature() (*ILS2T method*), 157
 get_timestamps() (*RTO1024 method*), 149
 get_ui_lower_limits() (*PSI9000 method*), 63
 get_uip_upper_limits() (*PSI9000 method*), 63
 get_voltage_current_setpoint() (*PSI9000 method*), 63
 GetAttr (class in *hvl_ccb.utils.conversion.utils*), 195
 go_home() (*NewportSMC100PP method*), 128
 go_to_configuration() (*NewportSMC100PP method*), 128
 GREEN_NOT_READY (*SafetyStatus attribute*), 55
 GREEN_READY (*SafetyStatus attribute*), 55

H

HARDWARE (*CryLasLaser.RepetitionRates attribute*), 40
 has_safeground (*TiePieOscilloscopeChannelConfig property*), 177
 HEAD (*CryLasLaser.AnswersStatus attribute*), 39
 Heinzinger (class in *hvl_ccb.dev.heinzinger.device*), 93
 HeinzingerConfig (class in *hvl_ccb.dev.heinzinger.device*), 95
 HeinzingerDeviceNotRecognizedError, 92
 HeinzingerError, 92
 HeinzingerSerialCommunication (class in *hvl_ccb.dev.heinzinger.base*), 90
 HeinzingerSerialCommunicationConfig (class in *hvl_ccb.dev.heinzinger.base*), 90
 HeinzingerSetValueError, 92
 HIGH (*LabJack.DIOStatus attribute*), 105
 high_resolution (*FuGProbusVSetRegisters property*), 86
 home_search_polling_interval (*NewportSMC100PPConfig attribute*), 132
 home_search_timeout (*NewportSMC100PPConfig attribute*), 132
 home_search_type (*NewportSMC100PPConfig attribute*), 132
 home_search_velocity (*NewportSMC100PPConfig attribute*), 132

HOME_STARTED (*NewportSMC100PPSerialCommunication.ControllerError attribute*), 134
 HomeSwitch (*NewportSMC100PPConfig.HomeSearch attribute*), 131
 HomeSwitch_and_Index (*NewportSMC100PPConfig.HomeSearch attribute*), 131
 HOMING (*NewportStates attribute*), 137
 HOMING_FROM_RS232 (*NewportSMC100PP.StateMessages attribute*), 125
 HOMING_FROM_SMC (*NewportSMC100PP.StateMessages attribute*), 125
 HOMING_TIMEOUT (*NewportSMC100PP.MotorErrors attribute*), 124
 host (*ModbusTcpCommunicationConfig attribute*), 20
 host (*OpcUaCommunicationConfig attribute*), 23
 host (*TcpCommunicationConfig attribute*), 29
 host (*TelnetCommunicationConfig attribute*), 31
 host (*VisaCommunicationConfig attribute*), 34
 HT (*NewportConfigCommands attribute*), 123
 HUNDRED_VOLT (*ACVoltageRange attribute*), 74
 HUNDRED_VOLT (*DCCVoltageRange attribute*), 74
 hvl_ccb
 module, 204
 hvl_ccb.comm
 module, 35
 hvl_ccb.comm.base
 module, 11
 hvl_ccb.comm.labjack_1jm
 module, 16
 hvl_ccb.comm.modbus_tcp
 module, 19
 hvl_ccb.comm.opc
 module, 21
 hvl_ccb.comm.serial
 module, 24
 hvl_ccb.comm.tcp
 module, 28
 hvl_ccb.comm.telnet
 module, 30
 hvl_ccb.comm.visa
 module, 32
 hvl_ccb.configuration
 module, 200
 hvl_ccb.dev
 module, 192
 hvl_ccb.dev.base
 module, 187
 hvl_ccb.dev.crylas
 module, 46
 hvl_ccb.dev.crylas.crylas
 module, 35
 hvl_ccb.dev.cube

```

    module, 61
hvl_ccb.dev.cube.alarms
    module, 46
hvl_ccb.dev.cube.base
    module, 47
hvl_ccb.dev.cube.constants
    module, 53
hvl_ccb.dev.cube.earthing_stick
    module, 55
hvl_ccb.dev.cube.errors
    module, 56
hvl_ccb.dev.cube.picube
    module, 57
hvl_ccb.dev.cube.support
    module, 60
hvl_ccb.dev.cube.switches
    module, 60
hvl_ccb.dev.ea_psi9000
    module, 67
hvl_ccb.dev.ea_psi9000.ea_psi9000
    module, 62
hvl_ccb.dev.fluke884x
    module, 76
hvl_ccb.dev.fluke884x.base
    module, 67
hvl_ccb.dev.fluke884x.constants
    module, 72
hvl_ccb.dev.fluke884x.ranges
    module, 73
hvl_ccb.dev.fug
    module, 89
hvl_ccb.dev.fug.fug
    module, 76
hvl_ccb.dev.heinzinger
    module, 97
hvl_ccb.dev.heinzinger.base
    module, 90
hvl_ccb.dev.heinzinger.constants
    module, 92
hvl_ccb.dev.heinzinger.device
    module, 93
hvl_ccb.dev.heinzinger.mixin
    module, 96
hvl_ccb.dev.highland_t560
    module, 102
hvl_ccb.dev.highland_t560.base
    module, 97
hvl_ccb.dev.highland_t560.channel
    module, 100
hvl_ccb.dev.highland_t560.device
    module, 100
hvl_ccb.dev.labjack
    module, 110
hvl_ccb.dev.labjack.labjack
    module, 103
hvl_ccb.dev.lauda
    module, 118
hvl_ccb.dev.lauda.lauda
    module, 111
hvl_ccb.dev.mbw973
    module, 122
hvl_ccb.dev.mbw973.mbw973
    module, 118
hvl_ccb.dev.newport
    module, 137
hvl_ccb.dev.newport.newport
    module, 123
hvl_ccb.dev.pfeiffer_tpg
    module, 144
hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg
    module, 138
hvl_ccb.dev.protocols
    module, 146
hvl_ccb.dev.protocols.sources
    module, 145
hvl_ccb.dev.rs_rto1024
    module, 154
hvl_ccb.dev.rs_rto1024.rs_rto1024
    module, 146
hvl_ccb.dev.se_ils2t
    module, 162
hvl_ccb.dev.se_ils2t.se_ils2t
    module, 155
hvl_ccb.dev.sst_luminox
    module, 167
hvl_ccb.dev.sst_luminox.sst_luminox
    module, 163
hvl_ccb.dev.technix
    module, 174
hvl_ccb.dev.technix.base
    module, 168
hvl_ccb.dev.technix.device
    module, 171
hvl_ccb.dev.tiepie
    module, 185
hvl_ccb.dev.tiepie.base
    module, 174
hvl_ccb.dev.tiepie.channel
    module, 176
hvl_ccb.dev.tiepie.device
    module, 179
hvl_ccb.dev.tiepie.generator
    module, 180
hvl_ccb.dev.tiepie.oscilloscope
    module, 182
hvl_ccb.dev.tiepie.utils
    module, 185
hvl_ccb.dev.utils
    module, 186

```

```

    module, 190
hvl_ccb.dev.visa
    module, 190
hvl_ccb.error
    module, 202
hvl_ccb.experiment_manager
    module, 202
hvl_ccb.utils
    module, 200
hvl_ccb.utils.conversion
    module, 196
hvl_ccb.utils.conversion.map_range
    module, 192
hvl_ccb.utils.conversion.sensor
    module, 193
hvl_ccb.utils.conversion.unit
    module, 194
hvl_ccb.utils.conversion.utils
    module, 195
hvl_ccb.utils.enum
    module, 196
hvl_ccb.utils.poller
    module, 198
hvl_ccb.utils.typing
    module, 199
hvl_ccb.utils.validation
    module, 199
hysteresis_compensation (New-
    portSMC100PPConfig attribute), 132
|_
ID (FuGProbusIVCommands attribute), 82
identification (Fluke8845a property), 68
Identification_error (PfeifferTPG.SensorStatus at-
    tribute), 138
identifier (LJMCommunicationConfig attribute), 18
identify_device() (FuG method), 77
identify_device() (Heinzinger method), 94
identify_sensors() (PfeifferTPG method), 139
IKR (PfeifferTPG.SensorTypes attribute), 139
IKR11 (PfeifferTPG.SensorTypes attribute), 139
IKR9 (PfeifferTPG.SensorTypes attribute), 139
ILS2T (class in hvl_ccb.dev.se_il2t.se_il2t), 155
ILS2T.ActionsPtp (class in
    hvl_ccb.dev.se_il2t.se_il2t), 155
ILS2T.Mode (class in hvl_ccb.dev.se_il2t.se_il2t), 155
ILS2T.Ref16Jog (class in hvl_ccb.dev.se_il2t.se_il2t),
    155
ILS2T.State (class in hvl_ccb.dev.se_il2t.se_il2t), 156
ILS2TConfig (class in hvl_ccb.dev.se_il2t.se_il2t), 159
ILS2TError, 160
ILS2TModbusTcpCommunication (class in
    hvl_ccb.dev.se_il2t.se_il2t), 160
ILS2TModbusTcpCommunicationConfig (class in
    hvl_ccb.dev.se_il2t.se_il2t), 160
ILS2TRegAddr (class in hvl_ccb.dev.se_il2t.se_il2t),
    161
ILS2TRegDatatype (class in
    hvl_ccb.dev.se_il2t.se_il2t), 161
IMMEDIATE (TriggerSource attribute), 73
IMMEDIATELY (FuGRampModes attribute), 86
IMPULSE_140KV (PowerSetup attribute), 55
IMR (PfeifferTPG.SensorTypes attribute), 139
INACTIVE (CryLasLaser.AnswersStatus attribute), 39
INACTIVE (DoorStatus attribute), 53
INACTIVE (SwitchStatus attribute), 61
inhibit (Technix property), 171
init_attenuation (CryLasAttenuatorConfig attribute),
    37
init_monitored_nodes() (OpcUaCommunication
    method), 21
init_shutter_status (CryLasLaserConfig attribute),
    43
initialize() (NewportSMC100PP method), 128
INITIALIZED (ExperimentStatus attribute), 203
INITIALIZING (ExperimentStatus attribute), 204
INITIALIZING (SafetyStatus attribute), 55
initiate_trigger() (Fluke8845a method), 68
INPUT (FuGProbusVRegisterGroups attribute), 85
INPUT (GateMode attribute), 98
input_range (TiePieOscilloscopeChannelConfig prop-
    erty), 177
INSTALL (AutoInstallMode attribute), 97
INT32 (ILS2TRegDatatype attribute), 162
INT_SYNTHESIZER (TriggerMode attribute), 100
interface_type (PSI9000VisaCommunicationConfig
    attribute), 66
interface_type (RTO1024VisaCommunicationConfig
    attribute), 154
interface_type (VisaCommunicationConfig attribute),
    34
internal (LabJack.CjcType attribute), 104
INTERNAL (LaudaProRp245eConfig.ExtControlModeEnum
    attribute), 115
IO_SCANNING (ILS2TRegAddr attribute), 161
IoScanningModeValueError, 162
is_configdataclass (AsyncCommunicationProtocol-
    Config attribute), 14
is_configdataclass (BaseCubeConfiguration at-
    tribute), 50
is_configdataclass (CryLasAttenuatorConfig at-
    tribute), 37
is_configdataclass (CryLasLaserConfig attribute),
    43
is_configdataclass (EmptyConfig attribute), 189, 201
is_configdataclass (Fluke8845aConfig attribute), 70
is_configdataclass (FuGConfig attribute), 78

```

is_configdataclass (*HeinzingerConfig attribute*), 96
is_configdataclass (*ILS2TConfig attribute*), 159
is_configdataclass (*LaudaProRp245eConfig attribute*), 116
is_configdataclass (*LJMCommunicationConfig attribute*), 18
is_configdataclass (*LuminoxConfig attribute*), 164
is_configdataclass (*MBW973Config attribute*), 120
is_configdataclass (*ModbusTcpCommunicationConfig attribute*), 20
is_configdataclass (*NewportSMC100PPConfig attribute*), 132
is_configdataclass (*OpcUaCommunicationConfig attribute*), 23
is_configdataclass (*PfeifferTPGConfig attribute*), 141
is_configdataclass (*T560Config attribute*), 102
is_configdataclass (*TcpCommunicationConfig attribute*), 29
is_configdataclass (*TechnixConfig attribute*), 173
is_configdataclass (*TiePieDeviceConfig attribute*), 175
is_configdataclass (*VisaCommunicationConfig attribute*), 34
is_data_ready_polling_interval_sec (*TiePieDeviceConfig attribute*), 175
is_done() (*MBW973 method*), 119
is_error() (*ExperimentManager method*), 203
is_finished() (*ExperimentManager method*), 203
is_in_range() (*ILS2TRegDatatype method*), 162
is_inactive (*CryLasLaser.LaserStatus property*), 39
is_measurement_data_ready() (*TiePieOscilloscope method*), 183
is_measurement_running() (*TiePieOscilloscope method*), 183
is_open (*Client property*), 21
is_open (*LJMCommunication property*), 16
is_open (*OpcUaCommunication property*), 22
is_open (*SerialCommunication property*), 25
is_open (*TelnetCommunication property*), 30
is_polling() (*Poller method*), 198
is_ready (*CryLasLaser.LaserStatus property*), 39
is_running() (*ExperimentManager method*), 203
is_started (*Technix property*), 171
is_triggered() (*TiePieOscilloscope method*), 183
is_valid_scale_range_reversed_str() (*PfeifferTPGConfig.Model method*), 141

J

J (*LabJack.ThermocoupleType attribute*), 106
jerk_time (*NewportSMC100PPConfig attribute*), 132
JOG (*ILS2T.Mode attribute*), 155
jog_run() (*ILS2T method*), 158
jog_stop() (*ILS2T method*), 158

JOGGING (*NewportStates attribute*), 137
JOGGING_FROM_DISABLE (*NewportSMC100PP.StateMessages attribute*), 125
JOGGING_FROM_READY (*NewportSMC100PP.StateMessages attribute*), 125
JOGN_FAST (*ILS2TRegAddr attribute*), 161
JOGN_SLOW (*ILS2TRegAddr attribute*), 161
JR (*NewportConfigCommands attribute*), 123

K

K (*LabJack.TemperatureUnit attribute*), 105
K (*LabJack.ThermocoupleType attribute*), 106
K (*Temperature attribute*), 195
KELVIN (*Temperature attribute*), 195
keys() (*AsyncCommunicationProtocolConfig class method*), 14
keys() (*BaseCubeConfiguration class method*), 50
keys() (*BaseCubeOpcUaCommunicationConfig class method*), 52
keys() (*CryLasAttenuatorConfig class method*), 37
keys() (*CryLasAttenuatorSerialCommunicationConfig class method*), 38
keys() (*CryLasLaserConfig class method*), 43
keys() (*CryLasLaserSerialCommunicationConfig class method*), 45
keys() (*EmptyConfig class method*), 189, 201
keys() (*Fluke8845aConfig class method*), 70
keys() (*Fluke8845aTelnetCommunicationConfig class method*), 71
keys() (*FuGConfig class method*), 78
keys() (*FuGSerialCommunicationConfig class method*), 88
keys() (*HeinzingerConfig class method*), 96
keys() (*HeinzingerSerialCommunicationConfig class method*), 91
keys() (*ILS2TConfig class method*), 159
keys() (*ILS2TModbusTcpCommunicationConfig class method*), 160
keys() (*LaudaProRp245eConfig class method*), 116
keys() (*LaudaProRp245eTcpCommunicationConfig class method*), 117
keys() (*LJMCommunicationConfig class method*), 18
keys() (*LuminoxConfig class method*), 164
keys() (*LuminoxSerialCommunicationConfig class method*), 167
keys() (*MBW973Config class method*), 120
keys() (*MBW973SerialCommunicationConfig class method*), 122
keys() (*ModbusTcpCommunicationConfig class method*), 20
keys() (*NewportSMC100PPConfig class method*), 132

<code>keys()</code> (<i>NewportSMC100PPSerialCommunicationConfig class method</i>), 136		<code>LaudaProRp245eCommand</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 113
<code>keys()</code> (<i>OpcUaCommunicationConfig class method</i>), 23		<code>LaudaProRp245eCommandError</code> , 115
<code>keys()</code> (<i>PfeifferTPGConfig class method</i>), 141		<code>LaudaProRp245eConfig</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 115
<code>keys()</code> (<i>PfeifferTPGSerialCommunicationConfig class method</i>), 143		<code>LaudaProRp245eConfig.ExtControlModeEnum</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 115
<code>keys()</code> (<i>PICubeConfiguration class method</i>), 58		<code>LaudaProRp245eConfig.OperationModeEnum</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 115
<code>keys()</code> (<i>PICubeOpcUaCommunicationConfig class method</i>), 59	class	<code>LaudaProRp245eTcpCommunication</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 116
<code>keys()</code> (<i>PSI9000Config class method</i>), 65		<code>LaudaProRp245eTcpCommunicationConfig</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 117
<code>keys()</code> (<i>PSI9000VisaCommunicationConfig class method</i>), 66	class	<code>LEM4000S</code> (class <i>in hvl_ccb.utils.conversion.sensor</i>), 193
<code>keys()</code> (<i>RTO1024Config class method</i>), 153		<code>LF</code> (<i>FuGTerminators attribute</i>), 89
<code>keys()</code> (<i>RTO1024VisaCommunicationConfig class method</i>), 154	class	<code>LFCR</code> (<i>FuGTerminators attribute</i>), 89
<code>keys()</code> (<i>SerialCommunicationConfig class method</i>), 26		<code>LINE_1</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>T560CommunicationConfig class method</i>), 99		<code>LINE_10</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>T560Config class method</i>), 102		<code>LINE_11</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>TcpCommunicationConfig class method</i>), 29		<code>LINE_12</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>TechnixConfig class method</i>), 173		<code>LINE_13</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>TechnixSerialCommunicationConfig class method</i>), 169	class	<code>LINE_14</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>TechnixTelnetCommunicationConfig class method</i>), 170	class	<code>LINE_15</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>TelnetCommunicationConfig class method</i>), 31		<code>LINE_2</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>TiePieDeviceConfig class method</i>), 175		<code>LINE_3</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>VisaCommunicationConfig class method</i>), 34		<code>LINE_4</code> (<i>MessageBoard attribute</i>), 54
<code>keys()</code> (<i>VisaDeviceConfig class method</i>), 192		<code>LINE_5</code> (<i>MessageBoard attribute</i>), 54
L		<code>LINE_6</code> (<i>MessageBoard attribute</i>), 54
<code>LabJack</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 103		<code>LINE_7</code> (<i>MessageBoard attribute</i>), 54
<code>LabJack.AInRange</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 104	in	<code>LINE_8</code> (<i>MessageBoard attribute</i>), 54
<code>LabJack.BitLimit</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 104	in	<code>LINE_9</code> (<i>MessageBoard attribute</i>), 54
<code>LabJack.CalMicroAmpere</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 104	in	<code>list_devices()</code> (<i>TiePieOscilloscope static method</i>), 183
<code>LabJack.CjcType</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 104	in	<code>list_directory()</code> (<i>RTO1024 method</i>), 149
<code>LabJack.ClockFrequency</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 104	in	<code>LJMCommunication</code> (class <i>in hvl_ccb.comm.labjack_ljm</i>), 16
<code>LabJack.DeviceType</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 105	in	<code>LJMCommunicationConfig</code> (class <i>in hvl_ccb.comm.labjack_ljm</i>), 17
<code>LabJack.DIOStatus</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 105	in	<code>LJMCommunicationConfig.ConnectionType</code> (class <i>in hvl_ccb.comm.labjack_ljm</i>), 17
<code>LabJack.TemperatureUnit</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 105	in	<code>LJMCommunicationConfig.DeviceType</code> (class <i>in hvl_ccb.comm.labjack_ljm</i>), 17
<code>LabJack.ThermocoupleType</code> (class <i>in hvl_ccb.dev.labjack.labjack</i>), 105	in	<code>LJMCommunicationError</code> , 18
<code>LabJackError</code> , 110		<code>lm34</code> (<i>LabJack.CjcType attribute</i>), 104
<code>LabJackIdentifierDIOError</code> , 110		<code>LMT70A</code> (class <i>in hvl_ccb.utils.conversion.sensor</i>), 193
<code>laser_off()</code> (<i>CryLasLaser method</i>), 40		<code>load_configuration()</code> (<i>RTO1024 method</i>), 149
<code>laser_on()</code> (<i>CryLasLaser method</i>), 40		<code>load_device_configuration()</code> (<i>T560 method</i>), 101
<code>LaudaProRp245e</code> (class <i>in hvl_ccb.dev.lauda.lauda</i>), 111		<code>local_display()</code> (<i>RTO1024 method</i>), 149
		<code>LOCKED</code> (<i>DoorStatus attribute</i>), 53
		<code>LOW</code> (<i>LabJack.DIOStatus attribute</i>), 105
		<code>LOWER_TEMP</code> (<i>LaudaProRp245eCommand attribute</i>), 114
		<code>lower_temp</code> (<i>LaudaProRp245eConfig attribute</i>), 116
		<code>Luminox</code> (class <i>in hvl_ccb.dev.sst_luminox.sst_luminox</i>), 163

LuminoxConfig	(class <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 164	<i>in</i>	MBW973SerialCommunication	(class <i>hvl_ccb.dev.mbw973.mbw973</i>), 121	<i>in</i>
LuminoxError	, 165		MBW973SerialCommunicationConfig	(class <i>hvl_ccb.dev.mbw973.mbw973</i>), 121	<i>in</i>
LuminoxMeasurementType	(class <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 165	<i>in</i>	measure()	(<i>Fluke8845a method</i>), 68	
LuminoxMeasurementTypeDict	(in module <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 165		measure()	(<i>PfeifferTPG method</i>), 140	
LuminoxMeasurementTypeError	, 165		measure_all()	(<i>PfeifferTPG method</i>), 140	
LuminoxMeasurementTypeValue	(in module <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 165		measure_current()	(<i>DeprecatedHeinzingerMixin static method</i>), 96	
LuminoxOutputMode	(class <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 166	<i>in</i>	measure_voltage()	(<i>DeprecatedHeinzingerMixin static method</i>), 96	
LuminoxOutputModeError	, 166		measure_voltage_current()	(<i>PSI9000 method</i>), 63	
LuminoxSerialCommunication	(class <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 166	<i>in</i>	measurement_function	(<i>Fluke8845a property</i>), 69	
LuminoxSerialCommunicationConfig	(class <i>hvl_ccb.dev.sst_luminox.sst_luminox</i>), 166	<i>in</i>	MeasurementFunction	(class <i>hvl_ccb.dev.fluke884x.constants</i>), 72	<i>in</i>
LUT (LMT70A attribute)	, 194		MEDIUM_FILTER	(<i>FilterRange attribute</i>), 75	
M			MessageBoard	(class in <i>hvl_ccb.dev.cube.constants</i>), 53	
MANUAL (SwitchOperatingStatus attribute)	, 61		micro_step_per_full_step_factor	(<i>NewportSMC100PPConfig attribute</i>), 132	
MapBitAsymRange	(class <i>hvl_ccb.utils.conversion.map_range</i>), 192	<i>in</i>	MILLIMETER_MERCURY	(<i>Pressure attribute</i>), 194	
MapBitSymRange	(class <i>hvl_ccb.utils.conversion.map_range</i>), 193	<i>in</i>	MINIMUM	(<i>LabJack.ClockFrequency attribute</i>), 104	
MapRanges (class in <i>hvl_ccb.utils.conversion.map_range</i>), 193			MMHG	(<i>Pressure attribute</i>), 194	
MARK (SerialCommunicationParity attribute)	, 27		ModbusTcpCommunication	(class <i>hvl_ccb.comm.modbus_tcp</i>), 19	<i>in</i>
max_current (FuG property)	, 77		ModbusTcpCommunicationConfig	(class <i>hvl_ccb.comm.modbus_tcp</i>), 20	<i>in</i>
max_current (Heinzinger property)	, 94		ModbusTcpConnectionFailedError	, 21	
max_current (Source property)	, 145		model (PfeifferTPGConfig attribute)	, 141	
max_current (Technix property)	, 171		module		
max_current (TechnixConfig attribute)	, 173		<i>hvl_ccb</i> , 204		
max_current_hardware (FuG property)	, 77		<i>hvl_ccb.comm</i> , 35		
max_current_hardware (Heinzinger property)	, 94		<i>hvl_ccb.comm.base</i> , 11		
max_pr_number (LaudaProRp245eConfig attribute)	, 116		<i>hvl_ccb.comm.labjack_ljm</i> , 16		
max_pump_level (LaudaProRp245eConfig attribute)	, 116		<i>hvl_ccb.comm.modbus_tcp</i> , 19		
max_timeout_retry_nr (OpcUaCommunicationConfig attribute)	, 23		<i>hvl_ccb.comm.opc</i> , 21		
max_voltage (FuG property)	, 77		<i>hvl_ccb.comm.serial</i> , 24		
max_voltage (Heinzinger property)	, 94		<i>hvl_ccb.comm.tcp</i> , 28		
max_voltage (Source property)	, 145		<i>hvl_ccb.comm.telnet</i> , 30		
max_voltage (Technix property)	, 172		<i>hvl_ccb.comm.visa</i> , 32		
max_voltage (TechnixConfig attribute)	, 173		<i>hvl_ccb.configuration</i> , 200		
max_voltage_hardware (FuG property)	, 77		<i>hvl_ccb.dev</i> , 192		
max_voltage_hardware (Heinzinger property)	, 94		<i>hvl_ccb.dev.base</i> , 187		
MAXIMUM (LabJack.ClockFrequency attribute)	, 104		<i>hvl_ccb.dev.crylas</i> , 46		
MBW973 (class in <i>hvl_ccb.dev.mbw973.mbw973</i>), 118			<i>hvl_ccb.dev.crylas.crylas</i> , 35		
MBW973Config (class in <i>hvl_ccb.dev.mbw973.mbw973</i>), 120			<i>hvl_ccb.dev.cube</i> , 61		
MBW973ControlRunningError	, 121		<i>hvl_ccb.dev.cube.alarms</i> , 46		
MBW973Error	, 121		<i>hvl_ccb.dev.cube.base</i> , 47		
MBW973PumpRunningError	, 121		<i>hvl_ccb.dev.cube.constants</i> , 53		
			<i>hvl_ccb.dev.cube.earthing_stick</i> , 55		
			<i>hvl_ccb.dev.cube.errors</i> , 56		
			<i>hvl_ccb.dev.cube.picube</i> , 57		
			<i>hvl_ccb.dev.cube.support</i> , 60		
			<i>hvl_ccb.dev.cube.switches</i> , 60		
			<i>hvl_ccb.dev.ea_psi9000</i> , 67		

hvl_ccb.dev.ea_psi9000.ea_psi9000, 62
hvl_ccb.dev.fluke884x, 76
hvl_ccb.dev.fluke884x.base, 67
hvl_ccb.dev.fluke884x.constants, 72
hvl_ccb.dev.fluke884x.ranges, 73
hvl_ccb.dev.fug, 89
hvl_ccb.dev.fug.fug, 76
hvl_ccb.dev.heinzinger, 97
hvl_ccb.dev.heinzinger.base, 90
hvl_ccb.dev.heinzinger.constants, 92
hvl_ccb.dev.heinzinger.device, 93
hvl_ccb.dev.heinzinger.mixin, 96
hvl_ccb.dev.highland_t560, 102
hvl_ccb.dev.highland_t560.base, 97
hvl_ccb.dev.highland_t560.channel, 100
hvl_ccb.dev.highland_t560.device, 100
hvl_ccb.dev.labjack, 110
hvl_ccb.dev.labjack.labjack, 103
hvl_ccb.dev.lauda, 118
hvl_ccb.dev.lauda.lauda, 111
hvl_ccb.dev.mbw973, 122
hvl_ccb.dev.mbw973.mbw973, 118
hvl_ccb.dev.newport, 137
hvl_ccb.dev.newport.newport, 123
hvl_ccb.dev.pfeiffer_tpg, 144
hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg, 138
hvl_ccb.dev.protocols, 146
hvl_ccb.dev.protocols.sources, 145
hvl_ccb.dev.rs_rto1024, 154
hvl_ccb.dev.rs_rto1024.rs_rto1024, 146
hvl_ccb.dev.se_ilst2t, 162
hvl_ccb.dev.se_ilst2t.se_ilst2t, 155
hvl_ccb.dev.sst_luminox, 167
hvl_ccb.dev.sst_luminox.sst_luminox, 163
hvl_ccb.dev.technix, 174
hvl_ccb.dev.technix.base, 168
hvl_ccb.dev.technix.device, 171
hvl_ccb.dev.tiepie, 185
hvl_ccb.dev.tiepie.base, 174
hvl_ccb.dev.tiepie.channel, 176
hvl_ccb.dev.tiepie.device, 179
hvl_ccb.dev.tiepie.generator, 180
hvl_ccb.dev.tiepie.oscilloscope, 182
hvl_ccb.dev.tiepie.utils, 185
hvl_ccb.dev.utils, 190
hvl_ccb.dev.visa, 190
hvl_ccb.error, 202
hvl_ccb.experiment_manager, 202
hvl_ccb.utils, 200
hvl_ccb.utils.conversion, 196
hvl_ccb.utils.conversion.map_range, 192
hvl_ccb.utils.conversion.sensor, 193
hvl_ccb.utils.conversion.unit, 194

hvl_ccb.utils.conversion.utils, 195
hvl_ccb.utils.enum, 196
hvl_ccb.utils.poller, 198
hvl_ccb.utils.typing, 199
hvl_ccb.utils.validation, 199
MONITOR_I (*FuGProbusVRegisterGroups* attribute), 85
MONITOR_V (*FuGProbusVRegisterGroups* attribute), 85
most_recent_error (*FuGProbusVConfigRegisters* property), 83
motion_distance_per_full_step (*NewportSMC100PPConfig* attribute), 132
motor_config (*NewportSMC100PPConfig* property), 132
move_to_absolute_position() (*NewportSMC100PP* method), 128
move_to_relative_position() (*NewportSMC100PP* method), 129
move_wait_sec (*NewportSMC100PPConfig* attribute), 132
MOVING (*NewportSMC100PP.StateMessages* attribute), 125
MOVING (*NewportStates* attribute), 137
MS_NOMINAL_CURRENT (*PSI9000* attribute), 62
MS_NOMINAL_VOLTAGE (*PSI9000* attribute), 62
MULTI_COMMANDS_MAX (*VisaCommunication* attribute), 32
MULTI_COMMANDS_SEPARATOR (*VisaCommunication* attribute), 32

N

n_channels (*TiePieOscilloscope* property), 183
n_max_try_get_device (*TiePieDeviceConfig* attribute), 175
name (*Fluke8845aConfig* attribute), 70
NameEnum (class in *hvl_ccb.utils.enum*), 196
NAMES (*SerialCommunicationParity* attribute), 27
names() (*RTO1024.TriggerModes* class method), 146
namespace_index (*BaseCubeConfiguration* attribute), 50
NATIVEONLY (*TiePieOscilloscopeAutoResolutionModes* attribute), 184
NED_END_OF_TURN (*NewportSMC100PP.MotorErrors* attribute), 124
NEG (*ILS2T.Ref16Jog* attribute), 156
NEG_FAST (*ILS2T.Ref16Jog* attribute), 156
NEGATIVE (*FuGPolarities* attribute), 81
NEGATIVE (*Polarity* attribute), 54
negative_software_limit (*NewportSMC100PPConfig* attribute), 132
NewportConfigCommands (class in *hvl_ccb.dev.newport.newport*), 123
NewportControllerError, 123
NewportError, 124
NewportMotorError, 124

NewportMotorPowerSupplyWasCutError	124		noise_level_measurement_channel_1 (BaseCube-Configuration attribute), 51
NewportSerialCommunicationError	137	in	noise_level_measurement_channel_2 (BaseCube-Configuration attribute), 51
NewportSMC100PP (class hvl_ccb.dev.newport.newport)	124	in	noise_level_measurement_channel_3 (BaseCube-Configuration attribute), 51
NewportSMC100PP.MotorErrors (class hvl_ccb.dev.newport.newport)	124	in	noise_level_measurement_channel_4 (BaseCube-Configuration attribute), 51
NewportSMC100PP.StateMessages (class hvl_ccb.dev.newport.newport)	124	in	NONE (ILS2T.Ref16Jog attribute), 156
NewportSMC100PPConfig (class hvl_ccb.dev.newport.newport)	131	in	NONE (LabJack.ThermocoupleType attribute), 106
NewportSMC100PPConfig.EspStageConfig (class in hvl_ccb.dev.newport.newport)	131	in	None (PfeifferTPG.SensorTypes attribute), 139
NewportSMC100PPConfig.HomeSearch (class hvl_ccb.dev.newport.newport)	131	in	NONE (SerialCommunicationParity attribute), 27
NewportSMC100PPConfig (class in hvl_ccb.dev.newport.newport)	131	in	NORMAL (RTO1024.TriggerModes attribute), 146
NewportSMC100PPSerialCommunication (class hvl_ccb.dev.newport.newport)	133	in	noSen (PfeifferTPG.SensorTypes attribute), 139
NewportSMC100PPSerialCommunication.Controller (class in hvl_ccb.dev.newport.newport)	133	in	noSENSOR (PfeifferTPG.SensorTypes attribute), 139
NewportSMC100PPSerialCommunication		in	nr_trials_activate (LuminoxConfig attribute), 164
NewportSMC100PPSerialCommunication.Controller		in	HvLCommunicationProtocol (class in hvl_ccb.comm.base), 14
NewportSMC100PPSerialCommunicationConfig (class in hvl_ccb.dev.newport.newport)	135		Number (in module hvl_ccb.utils.typing), 199
NewportStates (class in hvl_ccb.dev.newport.newport)	137		number_of_decimals (HeinzingerConfig attribute), 96
NewportUncertainPositionError	137		number_of_recordings (Heinzinger property), 94
NO (FuGDigitalVal attribute)	79		number_of_sensors (PfeifferTPG property), 140
NO_ERROR (NewportSMC100PPSerialCommunication.Controller.ERROR attribute)	134	O	
NO_REF (NewportStates attribute)	137		ODD (SerialCommunicationParity attribute), 27
NO_REF_ESP_STAGE_ERROR (New-attribute), portSMC100PP.StateMessages	125		OFF (AutoInstallMode attribute), 97
NO_REF_FROM_CONFIG (New-attribute), portSMC100PP.StateMessages	125		OFF (FuGDigitalVal attribute), 79
NO_REF_FROM_DISABLED (New-attribute), portSMC100PP.StateMessages	125		OFF (GateMode attribute), 98
NO_REF_FROM_HOMING (New-attribute), portSMC100PP.StateMessages	125		OFF (TriggerMode attribute), 100
NO_REF_FROM_JOGGING (New-attribute), portSMC100PP.StateMessages	125		offset (TiePieGeneratorConfig property), 180
NO_REF_FROM_MOVING (New-attribute), portSMC100PP.StateMessages	125		OH (NewportConfigCommands attribute), 123
NO_REF_FROM_READY (New-attribute), portSMC100PP.StateMessages	125		Ok (PfeifferTPG.SensorStatus attribute), 138
NO_REF_FROM_RESET (New-attribute), portSMC100PP.StateMessages	125		on (FuG property), 77
No_sensor (PfeifferTPG.SensorStatus attribute)	138		ON (FuGDigitalVal attribute), 79
NO_SOURCE (PowerSetup attribute)	55		on (FuGProbusVDIRegisters property), 84
NOISE (TiePieGeneratorSignalType attribute)	181		ON (ILS2T.State attribute), 156
			ONE (LabJack.AInRange attribute), 104
			ONE (RecordingsEnum attribute), 93
			ONE (SerialCommunicationStopbits attribute), 27
			ONE_AMPERE (ACCurrentRange attribute), 73
			ONE_AMPERE (DCCurrentRange attribute), 74
			ONE_HUNDRED_MICRO_AMPERE (DCCurrentRange attribute), 74
			ONE_HUNDRED_MILLI_AMPERE (ACCurrentRange attribute), 73
			ONE_HUNDRED_MILLI_AMPERE (DCCurrentRange attribute), 74
			ONE_HUNDRED_MILLI_SECOND (ApertureRange attribute), 74
			ONE_HUNDRED_MILLI_VOLT (ACVoltageRange attribute), 74
			ONE_HUNDRED_MILLI_VOLT (DCVoltageRange attribute), 75

ONE_HUNDRED_MILLION_OHM (*ResistanceRange attribute*), 75
ONE_HUNDRED_OHM (*ResistanceRange attribute*), 75
ONE_HUNDRED_THOUSAND_OHM (*ResistanceRange attribute*), 75
ONE_HUNDREDTH (*LabJack.AInRange attribute*), 104
ONE_MILLI_AMPERE (*DCCurrentRange attribute*), 74
ONE_MILLION_OHM (*ResistanceRange attribute*), 75
ONE_POINT_FIVE (*SerialCommunicationStopbits attribute*), 27
ONE_SECOND (*ApertureRange attribute*), 74
ONE_TENTH (*LabJack.AInRange attribute*), 104
ONE_THOUSAND_OHM (*ResistanceRange attribute*), 75
ONE_THOUSAND_VOLT (*DCVoltageRange attribute*), 75
ONE_VOLT (*ACVoltageRange attribute*), 74
ONE_VOLT (*DCVoltageRange attribute*), 75
ONLYUPWARDSOFFTOZERO (*FuGRampModes attribute*), 86
OPC_MAX_YEAR (*BaseCube attribute*), 47
OPC_MIN_YEAR (*BaseCube attribute*), 47
OpcUaCommunication (*class in hvl_ccb.comm.opc*), 21
OpcUaCommunicationConfig (*class in hvl_ccb.comm.opc*), 22
OpcUaCommunicationIOError, 24
OpcUaCommunicationTimeoutError, 24
OpcUaSubHandler (*class in hvl_ccb.comm.opc*), 24
OPEN (*DoorStatus attribute*), 53
OPEN (*SwitchOperation attribute*), 61
OPEN (*SwitchStatus attribute*), 61
open() (*CommunicationProtocol method*), 14
open() (*LaudaProRp245eTcpCommunication method*), 117
open() (*LJMCommunication method*), 16
open() (*ModbusTcpCommunication method*), 19
open() (*NullCommunicationProtocol method*), 15
open() (*OpcUaCommunication method*), 22
open() (*SerialCommunication method*), 25
open() (*Tcp method*), 28
open() (*TelnetCommunication method*), 30
open() (*VisaCommunication method*), 33
open_interlock (*Technix property*), 172
open_shutter() (*CryLasLaser method*), 41
open_timeout (*VisaCommunicationConfig attribute*), 34
OPENED (*CryLasLaser.AnswersShutter attribute*), 39
OPENED (*CryLasLaserShutterStatus attribute*), 46
operate (*BaseCube property*), 48
operate (*PICube property*), 57
OPERATION_MODE (*LaudaProRp245eCommand attribute*), 114
operation_mode (*LaudaProRp245eConfig attribute*), 116
optional_defaults() (*AsyncCommunicationProtocolConfig class method*), 14
optional_defaults() (*BaseCubeConfiguration class method*), 51
optional_defaults() (*BaseCubeOpcUaCommunicationConfig class method*), 52
optional_defaults() (*CryLasAttenuatorConfig class method*), 37
optional_defaults() (*CryLasAttenuatorSerialCommunicationConfig class method*), 38
optional_defaults() (*CryLasLaserConfig class method*), 43
optional_defaults() (*CryLasLaserSerialCommunicationConfig class method*), 45
optional_defaults() (*EmptyConfig class method*), 189, 201
optional_defaults() (*Fluke8845aConfig class method*), 70
optional_defaults()
 (*Fluke8845aTelnetCommunicationConfig class method*), 71
optional_defaults() (*FuGConfig class method*), 78
optional_defaults() (*FuGSerialCommunicationConfig class method*), 88
optional_defaults() (*HeinzingerConfig class method*), 96
optional_defaults() (*HeinzingerSerialCommunicationConfig class method*), 91
optional_defaults() (*ILS2TConfig class method*), 159
optional_defaults()
 (*ILS2TModbusTcpCommunicationConfig class method*), 161
optional_defaults() (*LaudaProRp245eConfig class method*), 116
optional_defaults() (*LaudaProRp245eTcpCommunicationConfig class method*), 117
optional_defaults() (*LJMCommunicationConfig class method*), 18
optional_defaults() (*LuminoxConfig class method*), 164
optional_defaults() (*LuminoxSerialCommunicationConfig class method*), 167
optional_defaults() (*MBW973Config class method*), 120
optional_defaults()
 (*MBW973SerialCommunicationConfig class method*), 122
optional_defaults() (*ModbusTcpCommunicationConfig class method*), 20
optional_defaults() (*NewportSMC100PPConfig class method*), 132
optional_defaults() (*NewportSMC100PPSerialCommunicationConfig class method*), 136

optional_defaults() (*OpcUaCommunicationConfig class method*), 23
optional_defaults() (*PfeifferTPGConfig class method*), 141
optional_defaults() (*PfeifferTPGSerialCommunicationConfig class method*), 143
optional_defaults() (*PICubeConfiguration class method*), 58
optional_defaults() (*PICubeOpcUaCommunicationConfig class method*), 59
optional_defaults() (*PSI9000Config class method*), 65
optional_defaults() (*PSI9000VisaCommunicationConfig class method*), 66
optional_defaults() (*RTO1024Config class method*), 153
optional_defaults() (*RTO1024VisaCommunicationConfig class method*), 154
optional_defaults() (*SerialCommunicationConfig class method*), 26
optional_defaults() (*T560CommunicationConfig class method*), 99
optional_defaults() (*T560Config class method*), 102
optional_defaults() (*TcpCommunicationConfig class method*), 29
optional_defaults() (*TechnixConfig class method*), 173
optional_defaults() (*TechnixSerialCommunicationConfig class method*), 169
optional_defaults() (*TechnixTelnetCommunicationConfig class method*), 170
optional_defaults() (*TelnetCommunicationConfig class method*), 31
optional_defaults() (*TiePieDeviceConfig class method*), 175
optional_defaults() (*VisaCommunicationConfig class method*), 34
optional_defaults() (*VisaDeviceConfig class method*), 192
OSCILLOSCOPE (*TiePieDeviceType attribute*), 175
OT (*NewportConfigCommands attribute*), 123
out (*FuGProbusVDORegisters property*), 84
OUTPUT (*FuGProbusIVCommands attribute*), 82
OUTPUT (*GateMode attribute*), 98
output (*Heinzinger property*), 95
output (*Source property*), 145
output (*Technix property*), 172
output_off() (*DeprecatedHeinzingerMixin static method*), 97
output_off() (*FuGProbusIV method*), 82
output_on() (*DeprecatedHeinzingerMixin static method*), 97
OUTPUT_POWER_EXCEEDED (*NewportSMC100PP.MotorErrors attribute*), 124
OUTPUTONCMD (*FuGProbusVRegisterGroups attribute*), 85
OUTPUTX0 (*FuGProbusVRegisterGroups attribute*), 85
OUTPUTX1 (*FuGProbusVRegisterGroups attribute*), 85
OUTPUTX2 (*FuGProbusVRegisterGroups attribute*), 85
OUTPUTXCMD (*FuGProbusVRegisterGroups attribute*), 85
outX0 (*FuG property*), 77
outX1 (*FuG property*), 77
outX2 (*FuG property*), 77
outXCMD (*FuG property*), 77
Overrange (*PfeifferTPG.SensorStatus attribute*), 138

P

PA (*Pressure attribute*), 194
PARAM_MISSING_OR_INVALID (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 134
parity (*CryLasAttenuatorSerialCommunicationConfig attribute*), 38
parity (*CryLasLaserSerialCommunicationConfig attribute*), 45
parity (*FuGSerialCommunicationConfig attribute*), 88
parity (*HeinzingerSerialCommunicationConfig attribute*), 91
parity (*LuminoxSerialCommunicationConfig attribute*), 167
parity (*MBW973SerialCommunicationConfig attribute*), 122
parity (*NewportSMC100PPSerialCommunicationConfig attribute*), 136
parity (*PfeifferTPGSerialCommunicationConfig attribute*), 143
Parity (*SerialCommunicationConfig attribute*), 26
parity (*SerialCommunicationConfig attribute*), 27
parse_read_measurement_value() (*LuminoxMeasurementType method*), 165
partial_pressure_o2 (*LuminoxMeasurementType attribute*), 165
PASCAL (*Pressure attribute*), 194
pause() (*LaudaProRp245e method*), 111
pause_ramp() (*LaudaProRp245e method*), 111
PBR (*PfeifferTPG.SensorTypes attribute*), 139
PEAK_CURRENT_LIMIT (*NewportSMC100PP.MotorErrors attribute*), 124
peak_output_current_limit (*NewportSMC100PPConfig attribute*), 132
percent_o2 (*LuminoxMeasurementType attribute*), 165
PERIOD (*MeasurementFunction attribute*), 72
period (*T560 property*), 101
period_aperture (*Fluke8845a attribute*), 69

PfeifferTPG	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 138	<i>in</i>	<i>portSMC100PPSerialCommunication.ControllerErrors</i> attribute), 134
PfeifferTPG.SensorStatus	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 138	<i>in</i>	POSITIVE (<i>FuGPolarities</i> attribute), 81
PfeifferTPG.SensorTypes	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 138	<i>in</i>	POSITIVE (<i>Polarity</i> attribute), 54
PfeifferTPGConfig	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 141	<i>in</i>	<i>positive_software_limit</i> (New- <i>portSMC100PPConfig</i> attribute), 133
PfeifferTPGConfig.Model	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 141	<i>in</i>	<i>post_force_value()</i> (Newport <i>SMC100PPConfig</i> method), 133
PfeifferTPGError	, 142	<i>in</i>	<i>post_stop_pause_sec</i> (<i>TechnixConfig</i> attribute), 173
PfeifferTPGSerialCommunication	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 142	<i>in</i>	POUNDS_PER_SQUARE_INCH (<i>Pressure</i> attribute), 194
PfeifferTPGSerialCommunicationConfig	(class <i>hvl_ccb.dev.pfeiffer_tpg.pfeiffer_tpg</i>), 142	<i>in</i>	POWER_INVERTER_220V (<i>PowerSetup</i> attribute), 55
PICube	(class in <i>hvl_ccb.dev.cube.picube</i>), 57	<i>in</i>	<i>power_limit</i> (<i>PSI9000Config</i> attribute), 65
PICubeConfiguration	(class <i>hvl_ccb.dev.cube.picube</i>), 58	<i>in</i>	<i>power_setup</i> (<i>PICube</i> property), 58
PICubeOpcUaCommunication	(class <i>hvl_ccb.dev.cube.picube</i>), 59	<i>in</i>	<i>PowerSetup</i> (class in <i>hvl_ccb.dev.cube.constants</i>), 54
PICubeOpcUaCommunicationConfig	(class <i>hvl_ccb.dev.cube.picube</i>), 59	<i>in</i>	<i>pre_sample_ratio</i> (<i>TiePieOscilloscopeConfig</i> prop- erty), 184
PICubeTestParameterError	, 56	<i>in</i>	<i>prepare_ultra_segmentation()</i> (RTO1024 method), 149
PKR (PfeifferTPG.SensorTypes attribute)	, 139	<i>in</i>	<i>preserve_type()</i> (in module <i>hvl_ccb.utils.conversion.utils</i>), 195
Polarity (class in <i>hvl_ccb.dev.cube.constants</i>)	, 54	<i>in</i>	<i>Pressure</i> (class in <i>hvl_ccb.utils.conversion.unit</i>), 194
Polarity (class in <i>hvl_ccb.dev.highland_t560.base</i>)	, 98	<i>in</i>	<i>probe_offset</i> (<i>TiePieOscilloscopeChannelConfig</i> prop- erty), 177
POLARITY (FuGProbusIVCommands attribute)	, 82	<i>in</i>	<i>PSI</i> (<i>Pressure</i> attribute), 195
polarity (PICube property)	, 58	<i>in</i>	<i>PSI9000</i> (class in <i>hvl_ccb.dev.ea_psi9000.ea_psi9000</i>), 62
Poller (class in <i>hvl_ccb.utils.poller</i>)	, 198	<i>in</i>	<i>PSI9000Config</i> (class <i>hvl_ccb.dev.ea_psi9000.ea_psi9000</i>), 64
polling (LuminoxOutputMode attribute)	, 166	<i>in</i>	<i>PSI9000Error</i> , 65
polling_delay_sec (BaseCubeConfiguration attribute)	, 51	<i>at-</i>	<i>PSI9000VisaCommunication</i> (class <i>hvl_ccb.dev.ea_psi9000.ea_psi9000</i>), 66
polling_interval (MBW973Config attribute)	, 120	<i>at-</i>	<i>PSI9000VisaCommunicationConfig</i> (class <i>hvl_ccb.dev.ea_psi9000.ea_psi9000</i>), 66
polling_interval_sec (BaseCubeConfiguration attribute)	, 51	<i>at-</i>	<i>PT100</i> (<i>LabJack.ThermocoupleType</i> attribute), 106
polling_interval_sec (TechnixConfig attribute)	, 173	<i>at-</i>	<i>PT1000</i> (<i>LabJack.ThermocoupleType</i> attribute), 106
polling_period (CryLasLaserConfig attribute)	, 43	<i>at-</i>	<i>PT500</i> (<i>LabJack.ThermocoupleType</i> attribute), 106
polling_timeout (CryLasLaserConfig attribute)	, 43	<i>at-</i>	<i>PTP</i> (ILS2T.Mode attribute), 155
port (Fluke8845aTelnetCommunicationConfig attribute)	, 71	<i>at-</i>	<i>PublicPropertiesReprMixin</i> (class <i>hvl_ccb.dev.tiepie.utils</i>), 185
port (ModbusTcpCommunicationConfig attribute)	, 20	<i>at-</i>	<i>PULSE</i> (<i>TiePieGeneratorSignalType</i> attribute), 181
port (OpcUaCommunicationConfig attribute)	, 23	<i>at-</i>	<i>pump_init</i> (<i>LaudaProRp245eConfig</i> attribute), 116
port (SerialCommunicationConfig attribute)	, 27	<i>at-</i>	<i>PUMP_LEVEL</i> (<i>LaudaProRp245eCommand</i> attribute), 114
port (T560CommunicationConfig attribute)	, 99		
port (TcpCommunicationConfig attribute)	, 29		
port (TechnixTelnetCommunicationConfig attribute)	, 170		
port (TelnetCommunicationConfig attribute)	, 31		
port (VisaCommunicationConfig attribute)	, 34		
POS (ILS2T.Ref16Jog attribute)	, 156		
POS_END_OF_TURN (Newport <i>SMC100PP.MotorErrors</i> attribute)	, 124		
POS_FAST (ILS2T.Ref16Jog attribute)	, 156		
POSITION (ILS2TRegAddr attribute)	, 161		
POSITION_OUT_OF_LIMIT		(New-	

Q

QIL (NewportConfigCommands attribute)	, 123
QUERY (FuGProbusIVCommands attribute)	, 82
query() (CryLasLaserSerialCommunication method)	, 44
query() (Fluke8845aTelnetCommunication method)	, 70
query() (FuGSerialCommunication method)	, 87
query() (Newport <i>SMC100PPSerialCommunication</i> method)	, 134
query() (PfeifferTPGSerialCommunication method)	, 142

query() (*SyncCommunicationProtocol method*), 15
query() (*T560Communication method*), 98
query() (*VisaCommunication method*), 33
query_all() (*CryLasLaserSerialCommunication method*), 44
query_command() (*LaudaProRp245eTcpCommunication method*), 117
query_multiple() (*NewportSMC100PPSerialCommunication method*), 134
query_polling() (*Luminox method*), 163
query_status() (*Technix method*), 172
QUEUE (*AutoInstallMode attribute*), 98
QUICK_STOP (*SafetyStatus attribute*), 55
QUICKSTOP (*ILS2T.State attribute*), 156
quickstop() (*ILS2T method*), 158
quit_error() (*BaseCube method*), 49

R

R (*LabJack.ThermocoupleType attribute*), 106
raise_() (*FuGErrorcodes method*), 80
RAMP_ACC (*ILS2TRegAddr attribute*), 161
RAMP_CONTINUE (*LaudaProRp245eCommand attribute*), 114
RAMP_DECEL (*ILS2TRegAddr attribute*), 161
RAMP_DELETE (*LaudaProRp245eCommand attribute*), 114
RAMP_ITERATIONS (*LaudaProRp245eCommand attribute*), 114
RAMP_N_MAX (*ILS2TRegAddr attribute*), 161
RAMP_PAUSE (*LaudaProRp245eCommand attribute*), 114
RAMP_SELECT (*LaudaProRp245eCommand attribute*), 114
RAMP_SET (*LaudaProRp245eCommand attribute*), 114
RAMP_START (*LaudaProRp245eCommand attribute*), 114
RAMP_STOP (*LaudaProRp245eCommand attribute*), 114
RAMP_TYPE (*ILS2TRegAddr attribute*), 161
rampmode (*FuGProbusVSetRegisters property*), 86
ramprate (*FuGProbusVSetRegisters property*), 86
rampstate (*FuGProbusVSetRegisters property*), 86
RAMPUPWARDS (*FuGRampModes attribute*), 86
RangeEnum (*class in hvl_ccb.utils.enum*), 197
RATEDCURRENT (*FuGReadbackChannels attribute*), 87
RATEDVOLTAGE (*FuGReadbackChannels attribute*), 87
read() (*AsyncCommunicationProtocol method*), 11
read() (*BaseCube method*), 49
read() (*CryLasLaserSerialCommunication method*), 44
read() (*LaudaProRp245eTcpCommunication method*), 117
read() (*MBW973 method*), 119
read() (*OpcUaCommunication method*), 22
read() (*Tcp method*), 28
read_all() (*AsyncCommunicationProtocol method*), 12

read_bytes() (*AsyncCommunicationProtocol method*), 12
read_bytes() (*SerialCommunication method*), 25
read_bytes() (*TelnetCommunication method*), 30
read_float() (*MBW973 method*), 119
read_holding_registers() (*ModbusTcpCommunication method*), 19
read_input_registers() (*ModbusTcpCommunication method*), 19
read_int() (*MBW973 method*), 119
read_measurement() (*RTO1024 method*), 149
read_measurements() (*MBW973 method*), 119
read_name() (*LJMCommunication method*), 16
read_nonempty() (*AsyncCommunicationProtocol method*), 12
read_output_while_polling (*TechnixConfig attribute*), 173
read_resistance() (*LabJack method*), 108
read_single_bytes() (*SerialCommunication method*), 25
read_streaming() (*Luminox method*), 164
read_termination (*VisaCommunicationConfig attribute*), 35
read_text() (*AsyncCommunicationProtocol method*), 12
read_text() (*NewportSMC100PPSerialCommunication method*), 135
read_text_nonempty() (*AsyncCommunicationProtocol method*), 12
READ_TEXT_SKIP_PREFIXES (*CryLasLaserSerialCommunication attribute*), 44
read_thermocouple() (*LabJack method*), 108
readback_data (*FuGProbusVConfigRegisters property*), 83
READBACKCHANNEL (*FuGProbusIVCommands attribute*), 82
ready (*BaseCube property*), 49
READY (*CryLasLaser.AnswersStatus attribute*), 39
READY (*ILS2T.State attribute*), 156
READY (*NewportStates attribute*), 137
READY_ACTIVE (*CryLasLaser.LaserStatus attribute*), 39
READY_FROM_DISABLE (*NewportSMC100PP.StateMessages attribute*), 125
READY_FROM_HOMING (*NewportSMC100PP.StateMessages attribute*), 125
READY_FROM_JOGGING (*NewportSMC100PP.StateMessages attribute*), 125
READY_FROM_MOVING (*NewportSMC100PP.StateMessages attribute*), 125
READY_INACTIVE (*CryLasLaser.LaserStatus attribute*),

39
record_length (*TiePieOscilloscopeConfig* property), 184
RecordingsEnum (class in *hvl_ccb.dev.heinzinger.constants*), 92
RED_OPERATE (*SafetyStatus* attribute), 55
RED_READY (*SafetyStatus* attribute), 55
reg_3 (*FuGProbusVDIRegisters* property), 84
RegAddr (*ILS2T* attribute), 156
RegDatatype (*ILS2T* attribute), 156
register_pulse_time (*TechnixConfig* attribute), 173
RELATIVE (*TiePieOscilloscopeTriggerLevelMode* attribute), 178
RELATIVE_POSITION_MOTOR (*ILS2T.ActionsPtp* attribute), 155
RELATIVE_POSITION_TARGET (*ILS2T.ActionsPtp* attribute), 155
remote (*Technix* property), 172
remove_device() (*DeviceSequenceMixin* method), 188
require_block_measurement_support (*TiePieDeviceConfig* attribute), 175
required_keys() (*AsyncCommunicationProtocolConfig* class method), 14
required_keys() (*BaseCubeConfiguration* class method), 51
required_keys() (*BaseCubeOpcUaCommunicationConfig* class method), 52
required_keys() (*CryLasAttenuatorConfig* class method), 37
required_keys() (*CryLasAttenuatorSerialCommunicationConfig* class method), 38
required_keys() (*CryLasLaserConfig* class method), 43
required_keys() (*CryLasLaserSerialCommunicationConfig* class method), 45
required_keys() (*EmptyConfig* class method), 189, 201
required_keys() (*Fluke8845aConfig* class method), 70
required_keys() (*Fluke8845aTelnetCommunicationConfig* class method), 71
required_keys() (*FuGConfig* class method), 78
required_keys() (*FuGSerialCommunicationConfig* class method), 88
required_keys() (*HeinzingerConfig* class method), 96
required_keys() (*HeinzingerSerialCommunicationConfig* class method), 91
required_keys() (*ILS2TConfig* class method), 160
required_keys() (*ILS2TModbusTcpCommunicationConfig* class method), 161
required_keys() (*LaudaProRp245eConfig* class method), 116
required_keys() (*LaudaProRp245eTcpCommunicationConfig* class method), 118
required_keys() (*LJMCommunicationConfig* class method), 18
required_keys() (*LuminoxConfig* class method), 164
required_keys() (*LuminoxSerialCommunicationConfig* class method), 167
required_keys() (*MBW973Config* class method), 121
required_keys() (*MBW973SerialCommunicationConfig* class method), 122
required_keys() (*ModbusTcpCommunicationConfig* class method), 20
required_keys() (*NewportSMC100PPConfig* class method), 133
required_keys() (*NewportSMC100PPSerialCommunicationConfig* class method), 136
required_keys() (*OpcUaCommunicationConfig* class method), 23
required_keys() (*PfeifferTPGConfig* class method), 142
required_keys() (*PfeifferTPGSerialCommunicationConfig* class method), 143
required_keys() (*PICubeConfiguration* class method), 59
required_keys() (*PICubeOpcUaCommunicationConfig* class method), 60
required_keys() (*PSI9000Config* class method), 65
required_keys() (*PSI9000VisaCommunicationConfig* class method), 66
required_keys() (*RTO1024Config* class method), 153
required_keys() (*RTO1024VisaCommunicationConfig* class method), 154
required_keys() (*SerialCommunicationConfig* class method), 27
required_keys() (*T560CommunicationConfig* class method), 99
required_keys() (*T560Config* class method), 102
required_keys() (*TcpCommunicationConfig* class method), 29
required_keys() (*TechnixConfig* class method), 173
required_keys() (*TechnixSerialCommunicationConfig* class method), 169
required_keys() (*TechnixTelnetCommunicationConfig* class method), 170
required_keys() (*TelnetCommunicationConfig* class method), 31
required_keys() (*TiePieDeviceConfig* class method), 175
required_keys() (*VisaCommunicationConfig* class method), 35
required_keys() (*VisaDeviceConfig* class method), 192
RESET (*FuGProbusIVCommands* attribute), 82
reset() (*Fluke8845a* method), 69
reset() (*FuGProbusIV* method), 82

reset() (*NewportSMC100PP method*), 129
reset() (*VisaDevice method*), 191
reset_error() (*ILS2T method*), 158
reset_interface() (*Heinzinger method*), 95
reset_ramp() (*LaudaProRp245e method*), 112
ResistanceRange (class) in *hvl_ccb.dev.fluke884x.ranges*, 75
resolution (*TiePieOscilloscopeConfig property*), 184
response_sleep_time (*CryLasAttenuatorConfig attribute*), 37
RISING (*TiePieOscilloscopeTriggerKind attribute*), 178
RMS_CURRENT_LIMIT (*NewportSMC100PP.MotorErrors attribute*), 124
rpm_max_init (*ILS2TConfig attribute*), 160
rs485_address (*NewportSMC100PPConfig attribute*), 133
RT01024 (class in *hvl_ccb.dev.rs_rto1024.rs_rto1024*), 146
RT01024.TriggerModes (class) in *hvl_ccb.dev.rs_rto1024.rs_rto1024*, 146
RT01024Config (class) in *hvl_ccb.dev.rs_rto1024.rs_rto1024*, 153
RT01024Error, 153
RT01024VisaCommunication (class) in *hvl_ccb.dev.rs_rto1024.rs_rto1024*, 153
RT01024VisaCommunicationConfig (class) in *hvl_ccb.dev.rs_rto1024.rs_rto1024*, 153
run() (*ExperimentManager method*), 203
run() (*LaudaProRp245e method*), 112
run_continuous_acquisition() (*RT01024 method*), 149
run_single_acquisition() (*RT01024 method*), 149
RUNNING (*ExperimentStatus attribute*), 204

S

S (*LabJack.ThermocoupleType attribute*), 106
SA (*NewportConfigCommands attribute*), 123
SafeGround (class in *hvl_ccb.dev.tiepie.channel*), 176
SafetyStatus (class in *hvl_ccb.dev.cube.constants*), 55
sample_frequency (*TiePieOscilloscopeConfig property*), 184
sample_rate (*TiePieOscilloscopeConfig property*), 184
save_configuration() (*RT01024 method*), 149
save_device_configuration() (*T560 method*), 101
save_waveform_history() (*RT01024 method*), 150
SCALE (*ILS2TRegAddr attribute*), 161
ScalingFactorValueError, 162
screw_scaling (*NewportSMC100PPConfig attribute*), 133
send_command() (*NewportSMC100PPSerialCommunication method*), 135
send_command() (*PfeifferTPGSerialCommunication method*), 142
send_hello() (*Client method*), 21
send_pulses() (*LabJack method*), 108
send_stop() (*NewportSMC100PPSerialCommunication method*), 135
Sensor (class in *hvl_ccb.utils.conversion.sensor*), 194
Sensor_error (*PfeifferTPG.SensorStatus attribute*), 138
Sensor_off (*PfeifferTPG.SensorStatus attribute*), 138
sensor_status (*LuminoxMeasurementType attribute*), 165
SERIAL (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 115
serial_number (*LuminoxMeasurementType attribute*), 165
serial_number (*TiePieDeviceConfig attribute*), 175
SerialCommunication (class in *hvl_ccb.comm.serial*), 24
SerialCommunicationBytesize (class) in *hvl_ccb.comm.serial*, 25
SerialCommunicationConfig (class) in *hvl_ccb.comm.serial*, 26
SerialCommunicationIOError, 27
SerialCommunicationParity (class) in *hvl_ccb.comm.serial*, 27
SerialCommunicationStopbits (class) in *hvl_ccb.comm.serial*, 27
Server (class in *hvl_ccb.comm.opc*), 24
set_acceleration() (*NewportSMC100PP method*), 129
set_acquire_length() (*RT01024 method*), 150
set_ain_differential() (*LabJack method*), 108
set_ain_range() (*LabJack method*), 108
set_ain_resolution() (*LabJack method*), 109
set_ain_thermocouple() (*LabJack method*), 109
set_analog_output() (*LabJack method*), 109
set_attenuation() (*CryLasAttenuator method*), 36
set_channel_offset() (*RT01024 method*), 150
set_channel_position() (*RT01024 method*), 150
set_channel_range() (*RT01024 method*), 151
set_channel_scale() (*RT01024 method*), 151
set_channel_state() (*RT01024 method*), 151
set_clock() (*LabJack method*), 109
set_control_mode() (*LaudaProRp245e method*), 112
set_current (*Heinzinger property*), 95
set_current (*Source property*), 145
set_current (*Technix property*), 172
set_digital_output() (*LabJack method*), 110
set_external_temp() (*LaudaProRp245e method*), 112
set_full_scale_mbar() (*PfeifferTPG method*), 140
set_full_scale_unitless() (*PfeifferTPG method*), 140
set_init_attenuation() (*CryLasAttenuator method*), 36

```

set_init_shutter_status() (CryLasLaser method), 41
set_jog_speed() (ILS2T method), 158
set_lower_limits() (PSI9000 method), 63
set_max_acceleration() (ILS2T method), 158
set_max_deceleration() (ILS2T method), 158
set_max_rpm() (ILS2T method), 158
set_measuring_options() (MBW973 method), 119
set_message_board() (BaseCube method), 49
set_motor_configuration() (NewportSMC100PP method), 129
set_negative_software_limit() (NewportSMC100PP method), 130
set_number_of_recordings() (Deprecated-HeinzingerMixin static method), 97
set_output() (PSI9000 method), 63
set_positive_software_limit() (NewportSMC100PP method), 130
set_pulse_energy() (CryLasLaser method), 41
set_pump_level() (LaudaProRp245e method), 112
set_ramp_iterations() (LaudaProRp245e method), 112
set_ramp_program() (LaudaProRp245e method), 112
set_ramp_segment() (LaudaProRp245e method), 113
set_ramp_type() (ILS2T method), 158
set_reference_point() (RTO1024 method), 151
set_register() (FuGProbusV method), 83
set_repetition_rate() (CryLasLaser method), 41
set_repetitions() (RTO1024 method), 152
set_status_board() (BaseCube method), 49
set_system_lock() (PSI9000 method), 64
set_temp_set_point() (LaudaProRp245e method), 113
set_transmission() (CryLasAttenuator method), 36
set_trigger_level() (RTO1024 method), 152
set_trigger_mode() (RTO1024 method), 152
set_trigger_source() (RTO1024 method), 152
set_upper_limits() (PSI9000 method), 64
set_voltage (Heinzinger property), 95
set_voltage (Source property), 145
set_voltage (Technix property), 172
set_voltage_current() (PSI9000 method), 64
SetAttr (class in hvl_ccb.utils.conversion.utils), 195
SETCURRENT (FuGProbusVRegisterGroups attribute), 85
setvalue (FuGProbusVSetRegisters property), 86
SETVOLTAGE (FuGProbusVRegisterGroups attribute), 85
SEVEN_HUNDRED_FIFTY_VOLT (ACVoltageRange attribute), 74
SEVENBITS (SerialCommunicationBytesize attribute), 26
SHORT_CIRCUIT (NewportSMC100PP.MotorErrors attribute), 124
shunt (LEM4000S attribute), 193
SHUTDOWN_CURRENT_LIMIT (PSI9000 attribute), 62
SHUTDOWN_VOLTAGE_LIMIT (PSI9000 attribute), 62
ShutterStatus (CryLasLaser attribute), 40
ShutterStatus (CryLasLaserConfig attribute), 42
signal_type (TiePieGeneratorConfig property), 180
SINE (TiePieGeneratorSignalType attribute), 181
SingleCommDevice (class in hvl_ccb.dev.base), 189
SIXBITS (SerialCommunicationBytesize attribute), 26
SIXTEEN (RecordingsEnum attribute), 93
SIXTEEN_BIT (TiePieOscilloscopeResolution attribute), 185
SL (NewportConfigCommands attribute), 123
SLOW_FILTER (FilterRange attribute), 75
SN (FuGReadbackChannels attribute), 87
SOFTWARE_INTERNAL_SIXTY (CryLasLaser.RepetitionRates attribute), 40
SOFTWARE_INTERNAL_TEN (CryLasLaser.RepetitionRates attribute), 40
SOFTWARE_INTERNAL_TWENTY (CryLasLaser.RepetitionRates attribute), 40
software_revision (LuminoxMeasurementType attribute), 165
Source (class in hvl_ccb.dev.protocols.sources), 145
SPACE (SerialCommunicationParity attribute), 27
SPECIALRAMPUPWARDS (FuGRampModes attribute), 86
spoll() (VisaCommunication method), 33
spoll_handler() (VisaDevice method), 191
SQUARE (TiePieGeneratorSignalType attribute), 181
SR (NewportConfigCommands attribute), 123
srq_mask (FuGProbusVConfigRegisters property), 83
srq_status (FuGProbusVConfigRegisters property), 83
stage_configuration (NewportSMC100PPConfig attribute), 133
START (LaudaProRp245eCommand attribute), 114
start() (BaseCube method), 49
start() (CryLasAttenuator method), 36
start() (CryLasLaser method), 41
start() (Device method), 187
start() (DeviceSequenceMixin method), 188
start() (ExperimentManager method), 203
start() (Fluke8845a method), 69
start() (FuG method), 78
start() (FuGProbusIV method), 82
start() (Heinzinger method), 95
start() (ILS2T method), 158
start() (LabJack method), 110
start() (LaudaProRp245e method), 113
start() (Luminox method), 164
start() (MBW973 method), 120
start() (NewportSMC100PP method), 130
start() (PfeifferTPG method), 140
start() (PSI9000 method), 64
start() (RTO1024 method), 152
start() (SingleCommDevice method), 190
start() (Technix method), 172
start() (TiePieGeneratorMixin method), 181

```

start() (*TiePieOscilloscope* method), 183
start() (*VisaDevice* method), 191
start_control() (*MBW973* method), 120
start_measurement() (*TiePieOscilloscope* method), 183
start_polling() (*Poller* method), 198
start_ramp() (*LaudaProRp245e* method), 113
STARTING (*ExperimentStatus* attribute), 204
States (*NewportSMC100PP* attribute), 125
status (*BaseCube* property), 49
status (*ExperimentManager* property), 203
status (*FuGProbusVConfigRegisters* property), 83
status (*FuGProbusVDORegisters* property), 85
status (*Technix* property), 172
STATUSBYTE (*FuGReadbackChannels* attribute), 87
STOP (*LaudaProRp245eCommand* attribute), 114
stop() (*BaseCube* method), 50
stop() (*CryLasLaser* method), 41
stop() (*Device* method), 187
stop() (*DeviceSequenceMixin* method), 188
stop() (*ExperimentManager* method), 203
stop() (*Fluke8845a* method), 69
stop() (*FuGProbusIV* method), 82
stop() (*Heinzinger* method), 95
stop() (*ILS2T* method), 158
stop() (*LabJack* method), 110
stop() (*LaudaProRp245e* method), 113
stop() (*Luminox* method), 164
stop() (*MBW973* method), 120
stop() (*NewportSMC100PP* method), 130
stop() (*PfeifferTPG* method), 141
stop() (*PSI9000* method), 64
stop() (*RTO1024* method), 152
stop() (*SingleCommDevice* method), 190
stop() (*Technix* method), 172
stop() (*TiePieGeneratorMixin* method), 181
stop() (*TiePieOscilloscope* method), 184
stop() (*VisaDevice* method), 191
stop_acquisition() (*RTO1024* method), 153
stop_measurement() (*TiePieOscilloscope* method), 184
stop_motion() (*NewportSMC100PP* method), 130
stop_polling() (*Poller* method), 198
stop_ramp() (*LaudaProRp245e* method), 113
STOP_SAFETY_STATUSES (in module *hvl_ccb.dev(cube.constants)*), 55
stopbits (*CryLasAttenuatorSerialCommunicationConfig* attribute), 38
stopbits (*CryLasLaserSerialCommunicationConfig* attribute), 45
stopbits (*FuGSerialCommunicationConfig* attribute), 88
stopbits (*HeinzingerSerialCommunicationConfig* attribute), 91
stopbits (*LuminoxSerialCommunicationConfig* attribute), 167
stopbits (*MBW973SerialCommunicationConfig* attribute), 122
stopbits (*NewportSMC100PPSerialCommunicationConfig* attribute), 137
stopbits (*PfeifferTPGSerialCommunicationConfig* attribute), 143
Stopbits (*SerialCommunicationConfig* attribute), 26
stopbits (*SerialCommunicationConfig* attribute), 27
streaming (*LuminoxOutputMode* attribute), 166
StrEnumBase (class in *hvl_ccb.utils.enum*), 197
sub_handler (*BaseCubeOpcUaCommunicationConfig* attribute), 52
sub_handler (*OpcUaCommunicationConfig* attribute), 23
SwitchOperatingStatus (class in *hvl_ccb.dev(cube.switches)*), 60
SwitchOperation (class in *hvl_ccb.dev(cube.switches)*), 61
SwitchOperationError, 56
SwitchStatus (class in *hvl_ccb.dev(cube.switches)*), 61
SyncCommunicationProtocol (class in *hvl_ccb.comm.base*), 15
SyncCommunicationProtocolConfig (class in *hvl_ccb.comm.base*), 15

T

T (*LabJack.ThermocoupleType* attribute), 106
t13_socket_1 (*BaseCube* attribute), 50
t13_socket_2 (*BaseCube* attribute), 50
t13_socket_3 (*BaseCube* attribute), 50
T1MS (*FuGMonitorModes* attribute), 80
T200MS (*FuGMonitorModes* attribute), 81
T20MS (*FuGMonitorModes* attribute), 81
T256US (*FuGMonitorModes* attribute), 81
T4 (*LabJack.DeviceType* attribute), 105
T4 (*LJMCommunicationConfig.DeviceType* attribute), 17
T40MS (*FuGMonitorModes* attribute), 81
T4MS (*FuGMonitorModes* attribute), 81
T560 (class in *hvl_ccb.dev.highland_t560.device*), 100
T560Communication (class in *hvl_ccb.dev.highland_t560.base*), 98
T560CommunicationConfig (class in *hvl_ccb.dev.highland_t560.base*), 98
T560Config (class in *hvl_ccb.dev.highland_t560.device*), 101
T560Error, 99
T7 (*LabJack.DeviceType* attribute), 105
T7 (*LJMCommunicationConfig.DeviceType* attribute), 17
T7_PRO (*LabJack.DeviceType* attribute), 105
T7_PRO (*LJMCommunicationConfig.DeviceType* attribute), 17
T800MS (*FuGMonitorModes* attribute), 81

T80MS (*FuGMonitorModes* attribute), 81
 target_pulse_energy (*CryLasLaser* property), 41
 Tcp (class in *hvl_ccb.comm.tcp*), 28
 TCP (*LJMCommunicationConfig.ConnectionType* attribute), 17
 TcpCommunicationConfig (class in *hvl_ccb.comm.tcp*), 28
 TCPIP_INSTR (*VisaCommunicationConfig.InterfaceType* attribute), 33
 TCPIP_SOCKET (*VisaCommunicationConfig.InterfaceType* attribute), 33
 TEC1 (*CryLasLaser.AnswersStatus* attribute), 39
 TEC2 (*CryLasLaser.AnswersStatus* attribute), 39
 Technix (class in *hvl_ccb.dev.technix.device*), 171
 TechnixConfig (class in *hvl_ccb.dev.technix.device*), 172
 TechnixError, 168
 TechnixFaultError, 168
 TechnixSerialCommunication (class in *hvl_ccb.dev.technix.base*), 168
 TechnixSerialCommunicationConfig (class in *hvl_ccb.dev.technix.base*), 168
 TechnixTelnetCommunication (class in *hvl_ccb.dev.technix.base*), 169
 TechnixTelnetCommunicationConfig (class in *hvl_ccb.dev.technix.base*), 169
 TelnetCommunication (class in *hvl_ccb.comm.telnet*), 30
 TelnetCommunicationConfig (class in *hvl_ccb.comm.telnet*), 31
 TelnetError, 32
 TEMP (*ILS2TRegAddr* attribute), 161
 TEMP_SET_POINT (*LaudaProRp245eCommand* attribute), 114
 temp_set_point_init (*LaudaProRp245eConfig* attribute), 116
 Temperature (class in *hvl_ccb.utils.conversion.unit*), 195
 temperature_sensor (*LuminoxMeasurementType* attribute), 165
 temperature_unit (*LMT70A* attribute), 194
 TEN (*LabJack.AInRange* attribute), 104
 TEN (*LabJack.CalMicroAmpere* attribute), 104
 TEN_AMPERE (*ACCurrrentRange* attribute), 73
 TEN_AMPERE (*DCCurrentRange* attribute), 74
 TEN_MHZ (*LabJack.ClockFrequency* attribute), 104
 TEN_MILLI_AMPERE (*ACCurrrentRange* attribute), 73
 TEN_MILLI_AMPERE (*DCCurrentRange* attribute), 74
 TEN_MILLI_SECOND (*ApertureRange* attribute), 74
 TEN_MILLION_OHM (*ResistanceRange* attribute), 75
 TEN_THOUSAND_OHM (*ResistanceRange* attribute), 75
 TEN_VOLT (*ACVoltageRange* attribute), 74
 TEN_VOLT (*DCVoltageRange* attribute), 75
 terminator (*AsyncCommunicationProtocolConfig* attribute), 14
 terminator (*CryLasAttenuatorSerialCommunicationConfig* attribute), 39
 terminator (*CryLasLaserSerialCommunicationConfig* attribute), 46
 terminator (*Fluke8845aTelnetCommunicationConfig* attribute), 71
 TERMINATOR (*FuGProbusIVCommands* attribute), 82
 terminator (*FuGProbusVConfigRegisters* property), 83
 terminator (*FuGSerialCommunicationConfig* attribute), 88
 terminator (*HeinzingerSerialCommunicationConfig* attribute), 92
 terminator (*LaudaProRp245eTcpCommunicationConfig* attribute), 118
 terminator (*LuminoxSerialCommunicationConfig* attribute), 167
 terminator (*MBW973SerialCommunicationConfig* attribute), 122
 terminator (*NewportSMC100PPSerialCommunicationConfig* attribute), 137
 terminator (*PfeifferTPGSerialCommunicationConfig* attribute), 144
 terminator (*T560CommunicationConfig* attribute), 99
 terminator_str() (*SerialCommunicationConfig* method), 27
 THIRTY_TWO_BIT (*LabJack.BitLimit* attribute), 104
 THREE_AMPERE (*ACCurrrentRange* attribute), 73
 THREE_AMPERE (*DCCurrentRange* attribute), 74
 TiePieDeviceConfig (class in *hvl_ccb.dev.tiepie.base*), 174
 TiePieDeviceType (class in *hvl_ccb.dev.tiepie.base*), 175
 TiePieError, 175
 TiePieGeneratorConfig (class in *hvl_ccb.dev.tiepie.generator*), 180
 TiePieGeneratorConfigLimits (class in *hvl_ccb.dev.tiepie.generator*), 180
 TiePieGeneratorMixin (class in *hvl_ccb.dev.tiepie.generator*), 180
 TiePieGeneratorSignalType (class in *hvl_ccb.dev.tiepie.generator*), 181
 TiePieHS5 (class in *hvl_ccb.dev.tiepie.device*), 179
 TiePieHS6 (class in *hvl_ccb.dev.tiepie.device*), 179
 TiePieOscilloscope (class in *hvl_ccb.dev.tiepie.oscilloscope*), 182
 TiePieOscilloscopeAutoResolutionModes (class in *hvl_ccb.dev.tiepie.oscilloscope*), 184
 TiePieOscilloscopeChannelConfig (class in *hvl_ccb.dev.tiepie.channel*), 176
 TiePieOscilloscopeChannelConfigLimits (class in *hvl_ccb.dev.tiepie.channel*), 177
 TiePieOscilloscopeChannelCoupling (class in

hvl_ccb.dev.tiepie.channel), 177
TiePieOscilloscopeConfig (class
hvl_ccb.dev.tiepie.oscilloscope), 184
TiePieOscilloscopeConfigLimits (class
hvl_ccb.dev.tiepie.oscilloscope), 185
TiePieOscilloscopeRange (class
hvl_ccb.dev.tiepie.channel), 178
TiePieOscilloscopeResolution (class
hvl_ccb.dev.tiepie.oscilloscope), 185
TiePieOscilloscopeTriggerKind (class
hvl_ccb.dev.tiepie.channel), 178
TiePieOscilloscopeTriggerLevelMode (class
hvl_ccb.dev.tiepie.channel), 178
TiePieWS5 (class in *hvl_ccb.dev.tiepie.device*), 179
timeout (*CryLasAttenuatorSerialCommunicationConfig*
attribute), 39
timeout (*CryLasLaserSerialCommunicationConfig*
attribute), 46
timeout (*FuGSerialCommunicationConfig* attribute), 88
timeout (*HeinzingerSerialCommunicationConfig*
attribute), 92
timeout (*LuminoxSerialCommunicationConfig* attribute), 167
timeout (*MBW973SerialCommunicationConfig* attribute), 122
timeout (*NewportSMC100PPSerialCommunicationConfig*
attribute), 137
timeout (*PfeifferTPGSerialCommunicationConfig*
attribute), 144
timeout (*SerialCommunicationConfig* attribute), 27
timeout (*TelnetCommunicationConfig* attribute), 31
timeout (*VisaCommunicationConfig* attribute), 35
timeout_interval (*BaseCubeConfiguration* attribute),
51
timeout_status_change (*BaseCubeConfiguration*
attribute), 51
timeout_test_parameters (*PICubeConfiguration*
attribute), 59
TORR (*Pressure* attribute), 195
TPG25xA (*PfeifferTPGConfig.Model* attribute), 141
TPGx6x (*PfeifferTPGConfig.Model* attribute), 141
TPR (*PfeifferTPG.SensorTypes* attribute), 139
transmission (*CryLasAttenuator* property), 36
TRIANGLE (*TiePieGeneratorSignalType* attribute), 181
trigger() (*Fluke8845a* method), 69
trigger_delay (*Fluke8845a* property), 69
trigger_enabled (*TiePieOscilloscopeChannelConfig*
property), 177
trigger_hysteresis (*TiePieOscilloscopeChannel*
Config property), 177
trigger_kind (*TiePieOscilloscopeChannelConfig*
property), 177
trigger_level (*T560* property), 101
trigger_level (*TiePieOscilloscopeChannelConfig*
property), 177
in **trigger_level_mode** (*TiePieOscilloscopeChannel*
Config property), 177
in **trigger_mode** (*T560* property), 101
trigger_source (*Fluke8845a* property), 69
in **trigger_timeout** (*TiePieOscilloscopeConfig* prop
erty), 185
in **TriggerMode** (class in *hvl_ccb.dev.highland_t560.base*),
99
in **TriggerSource** (class
hvl_ccb.dev.fluke884x.constants), 72
in **TWELVE_BIT** (*TiePieOscilloscopeResolution* attribute),
185
TWELVE_HUNDRED_FIFTY_KHZ (Lab
Jack.ClockFrequency attribute), 105
TWENTY_FIVE_HUNDRED_KHZ (LabJack.ClockFrequency
attribute), 105
TWENTY_MHZ (LabJack.ClockFrequency attribute), 105
TWENTY_VOLT (*TiePieOscilloscopeRange* attribute), 178
TWO (*RecordingsEnum* attribute), 93
TWO (*SerialCommunicationStopbits* attribute), 28
TWO_HUNDRED (LabJack.CalMicroAmpere attribute), 104
TWO_HUNDRED_MILLI_VOLT (*TiePieOscilloscopeRange*
attribute), 178
TWO_VOLT (*TiePieOscilloscopeRange* attribute), 178
TWO_WIRE_RESISTANCE (*MeasurementFunction* attribute), 72
two_wire_resistance_range (*Fluke8845a* attribute),
69

U

Underrange (*PfeifferTPG.SensorStatus* attribute), 138
Unit (class in *hvl_ccb.utils.conversion.unit*), 195
unit (*ILS2TModbusTcpCommunicationConfig* attribute),
161
unit (*ModbusTcpCommunicationConfig* attribute), 20
unit (*PfeifferTPG* property), 141
unit() (*RangeEnum* class method), 197
UNKNOWN (*TiePieGeneratorSignalType* attribute), 181
UNKNOWN (*TiePieOscilloscopeAutoResolutionModes*
attribute), 184
UNKNOWN (*TiePieOscilloscopeTriggerLevelMode*
attribute), 178
UNREADY_INACTIVE (*CryLasLaser.LaserStatus* attribute), 39
update_laser_status() (*CryLasLaser* method), 42
update_parameter (*OpcUaCommunicationConfig* attribute), 23
update_repetition_rate() (*CryLasLaser* method),
42
update_shutter_status() (*CryLasLaser* method), 42
update_target_pulse_energy() (*CryLasLaser*
method), 42

UpdateEspStageInfo (New-
portSMC100PPConfig.EspStageConfig attribute), 131

UPPER_TEMP (LaudaProRp245eCommand attribute), 114

upper_temp (LaudaProRp245eConfig attribute), 116

USB (LaudaProRp245eConfig.ExtControlModeEnum attribute), 115

USB (LJMCommunicationConfig.ConnectionType attribute), 17

use_external_clock() (T560 method), 101

user_position_offset (NewportSMC100PPConfig attribute), 133

user_steps() (ILS2T method), 159

V

VA (NewportConfigCommands attribute), 123

validate_and_resolve_host() (in module hvl_ccb.utils.validation), 199

validate_bool() (in module hvl_ccb.utils.validation), 199

validate_number() (in module hvl_ccb.utils.validation), 199

validate_pump_level() (LaudaProRp245e method), 113

validate_tcp_port() (in module hvl_ccb.utils.validation), 199

value (FuGProbusVMonitorRegisters property), 85

value (LabJack.AInRange property), 104

value_raw (FuGProbusVMonitorRegisters property), 85

ValueEnum (class in hvl_ccb.utils.enum), 197

VB (NewportConfigCommands attribute), 123

velocity (NewportSMC100PPConfig attribute), 133

visa_backend (VisaCommunicationConfig attribute), 35

VisaCommunication (class in hvl_ccb.comm.visa), 32

VisaCommunicationConfig (class in hvl_ccb.comm.visa), 33

VisaCommunicationConfig.InterfaceType (class in hvl_ccb.comm.visa), 33

VisaCommunicationError, 35

VisaDevice (class in hvl_ccb.dev.visa), 190

VisaDeviceConfig (class in hvl_ccb.dev.visa), 191

VOLT (ILS2TRegAddr attribute), 161

voltage (FuG property), 78

VOLTAGE (FuGProbusIVCommands attribute), 82

VOLTAGE (FuGReadbackChannels attribute), 87

voltage (Heinzinger property), 95

voltage (Source property), 145

voltage (Technix property), 172

VOLTAGE_AC (MeasurementFunction attribute), 72

voltage_actual (PICube property), 58

VOLTAGE_DC (MeasurementFunction attribute), 72

voltage_filter (Fluke8845a attribute), 69

voltage_lower_limit (PSI9000Config attribute), 65

voltage_max (PICube property), 58

voltage_monitor (FuG property), 78

voltage_primary (PICube property), 58

voltage_regulation (Technix property), 172

voltage_upper_limit (PSI9000Config attribute), 65

W

WAIT_AFTER_WRITE (VisaCommunication attribute), 32

wait_for_polling_result() (Poller method), 198

wait_operation_complete() (VisaDevice method), 191

wait_sec_initialisation (PSI9000Config attribute), 65

wait_sec_max_disable (ILS2TConfig attribute), 160

wait_sec_post_absolute_position (ILS2TConfig attribute), 160

wait_sec_post_activate (LuminoxConfig attribute), 165

wait_sec_post_cannot_disable (ILS2TConfig attribute), 160

wait_sec_post_enable (ILS2TConfig attribute), 160

wait_sec_post_relative_step (ILS2TConfig attribute), 160

wait_sec_pre_read_or_write (LaudaProRp245eTcpCommunicationConfig attribute), 118

wait_sec_read_text_nonempty (AsyncCommunicationProtocolConfig attribute), 14

wait_sec_read_text_nonempty (FuGSerialCommunicationConfig attribute), 88

wait_sec_read_text_nonempty (HeinzingerSerialCommunicationConfig attribute), 92

wait_sec_retry_get_device (TiePieDeviceConfig attribute), 175

wait_sec_settings_effect (PSI9000Config attribute), 65

wait_sec_stop_commands (FuGConfig attribute), 78

wait_sec_stop_commands (HeinzingerConfig attribute), 96

wait_sec_system_lock (PSI9000Config attribute), 65

wait_sec_trials_activate (LuminoxConfig attribute), 165

wait_timeout_retry_sec (OpcUaCommunicationConfig attribute), 23

wait_until_motor_initialized() (NewportSMC100PP method), 130

wait_until_ready() (CryLasLaser method), 42

waveform (TiePieGeneratorConfig property), 180

WIFI (LJMCommunicationConfig.ConnectionType attribute), 17

wrap.libtiepie_exception() (in module hvl_ccb.dev.tiepie.base), 176

write() (AsyncCommunicationProtocol method), 13

write() (BaseCube method), 50

write() (MBW973 method), 120

`write()` (*OpcUaCommunication method*), 22
`write()` (*Tcp method*), 28
`write()` (*VisaCommunication method*), 33
`write_absolute_position()` (*ILS2T method*), 159
`write_bytes()` (*AsyncCommunicationProtocol method*), 13
`write_bytes()` (*SerialCommunication method*), 25
`write_bytes()` (*TelnetCommunication method*), 30
`write_command()` (*Lau-daProRp245eTcpCommunication method*), 117
`write_name()` (*LJMCommunication method*), 17
`write_names()` (*LJMCommunication method*), 17
`write_registers()` (*ModbusTcpCommunication method*), 20
`write_relative_step()` (*ILS2T method*), 159
`write_termination` (*VisaCommunicationConfig attribute*), 35
`write_text()` (*AsyncCommunicationProtocol method*), 13
`WRONG_ESP_STAGE` (*NewportSMC100PP.MotorErrors attribute*), 124

X

`x_stat` (*FuGProbusVDIRegisters property*), 84
`XOUTPUTS` (*FuGProbusIVCommands attribute*), 82

Y

`YES` (*FuGDigitalVal attribute*), 79

Z

`ZX` (*NewportConfigCommands attribute*), 123