

---

# HVL Common Code Base Documentation

*Release 0.12.1*

**Mikolaj Rybiński, David Graber, Henrik Menne, Alise Chachereau,**

**Oct 31, 2022**



## CONTENTS:

<b>1</b>	<b>HVL Common Code Base</b>	<b>1</b>
1.1	Features . . . . .	1
1.2	Documentation . . . . .	5
1.3	Credits . . . . .	5
<b>2</b>	<b>Installation</b>	<b>7</b>
2.1	Stable release . . . . .	7
2.2	From sources . . . . .	7
2.3	Additional system libraries . . . . .	8
<b>3</b>	<b>Usage</b>	<b>9</b>
<b>4</b>	<b>API Documentation</b>	<b>11</b>
4.1	hvl_ccb . . . . .	11
<b>5</b>	<b>Contributing</b>	<b>187</b>
5.1	Types of Contributions . . . . .	187
5.2	Get Started! . . . . .	188
5.3	Merge Request Guidelines . . . . .	189
5.4	Tips . . . . .	189
5.5	Deploying . . . . .	190
<b>6</b>	<b>Credits</b>	<b>193</b>
6.1	Maintainers . . . . .	193
6.2	Authors . . . . .	193
6.3	Contributors . . . . .	193
<b>7</b>	<b>History</b>	<b>195</b>
7.1	0.12.1 (2022-10-31) . . . . .	195
7.2	0.12.0 (2022-10-17) . . . . .	195
7.3	0.11.1 (2022-09-15) . . . . .	195
7.4	0.11.0 (2022-06-22) . . . . .	196
7.5	0.10.3 (2022-03-21) . . . . .	196
7.6	0.10.2 (2022-02-28) . . . . .	196
7.7	0.10.1 (2022-01-24) . . . . .	196
7.8	0.10.0 (2022-01-17) . . . . .	197
7.9	0.9.0 (2022-01-07) . . . . .	197
7.10	0.8.5 (2021-11-05) . . . . .	198
7.11	0.8.4 (2021-10-22) . . . . .	198
7.12	0.8.3 (2021-09-27) . . . . .	198
7.13	0.8.2 (2021-08-27) . . . . .	198

7.14	0.8.1 (2021-08-13)	199
7.15	0.8.0 (2021-07-02)	199
7.16	0.7.1 (2021-06-04)	199
7.17	0.7.0 (2021-05-25)	199
7.18	0.6.1 (2021-05-08)	200
7.19	0.6.0 (2021-04-23)	200
7.20	0.5.0 (2020-11-11)	201
7.21	0.4.0 (2020-07-16)	201
7.22	0.3.5 (2020-02-18)	202
7.23	0.3.4 (2019-12-20)	202
7.24	0.3.3 (2019-05-08)	202
7.25	0.3.2 (2019-05-08)	202
7.26	0.3.1 (2019-05-02)	203
7.27	0.3 (2019-05-02)	203
7.28	0.2.1 (2019-04-01)	203
7.29	0.2.0 (2019-03-31)	203
7.30	0.1.0 (2019-02-06)	203

<b>8</b>	<b>Indices and tables</b>	<b>205</b>
	<b>Python Module Index</b>	<b>207</b>
	<b>Index</b>	<b>209</b>

---

**CHAPTER  
ONE**

---

## **HVL COMMON CODE BASE**

Python common code base (CCB) to control devices, which are used in high-voltage research. All implemented devices are used and tested in the High Voltage Laboratory ([HVL](#)) of the Federal Institute of Technology Zurich (ETH Zurich).

- Free software: GNU General Public License v3
- Copyright (c) 2019-2022 ETH Zurich, SIS ID and HVL D-ITET

### **1.1 Features**

For managing multi-device experiments instantiate the `ExperimentManager` utility class.

#### **1.1.1 Devices**

The device wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, various settings and options, as well as start/stop methods. Currently wrappers are available to control the following devices:

Function/Type	Devices
Bench Multimeter	Fluke 8845A and 8846A 6.5 Digit Precision Multimeter
Data acquisition	LabJack (T4, T7, T7-PRO; requires <a href="#">LJM Library</a> ) Pico Technology PT-104 Platinum Resistance Data Logger (requires <a href="#">PicoSDK/libusbpt104</a> )
Digital Delay Generator	Highland T560
Digital IO	LabJack (T4, T7, T7-PRO; requires <a href="#">LJM Library</a> )
Experiment control	HVL Cube with and without Power Inverter
Gas Analyser	MBW 973-SF6 gas dew point mirror analyzer Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges SST Luminox oxygen sensor
I2C host	TiePie (HS5, WS5; requires <a href="#">LibTiePie SDK</a> )
Laser	CryLaS pulsed laser CryLaS laser attenuator
Oscilloscope	Rhode & Schwarz RTO 1024 TiePie (HS5, HS6, WS5; requires <a href="#">LibTiePie SDK</a> )
Power supply	Elektro-Automatik PSI9000 FuG Elektronik Heinzinger PNC Technix capacitor charger
Stepper motor drive	Newport SMC100PP Schneider Electric ILS2T
Temperature control	
<b>2</b>	<b>Chapter 1. HVL Common Code Base</b> Lauda PRO RP 245 E circulation thermostat
Waveform generator	

Each device uses at least one standardised communication protocol wrapper.

### 1.1.2 Communication protocols

In hvl\_ccb by “communication protocol” we mean different levels of communication standards, from the low level actual communication protocols like serial communication to application level interfaces like VISA TCP standard. There are also devices in hvl\_ccb that use a dummy communication protocol; this is because these devices are build on proprietary manufacturer libraries that communicate with the corresponding devices, as in the case of TiePie or LabJack devices.

The communication protocol wrappers in hvl\_ccb provide a standardised API with configuration dataclasses, as well as open/close and read/write/query methods. Currently, wrappers for the following communication protocols are available:

Communication protocol	Devices using
Modbus TCP	Schneider Electric ILS2T stepper motor drive
OPC UA	HVL Cube with and without Power Inverter
Serial	CryLaS pulsed laser and laser attenuator FuG Elektronik power supply (e.g. capacitor charger HCK) using the Probus V protocol Heinzinger PNC power supply using Heinzinger Digital Interface I/II SST Luminox oxygen sensor MBW 973-SF6 gas dew point mirror analyzer Newport SMC100PP single axis driver for 2-phase stepper motors Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges Technix capacitor charger
TCP	Lauda PRO RP 245 E circulation thermostat
Telnet	Technix capacitor charger Fluke 8845A and 8846
VISA TCP	Elektro-Automatik PSI9000 DC power supply Rhode & Schwarz RTO 1024 oscilloscope
<i>proprietary</i>	LabJack (T4, T7, T7-PRO) devices, which communicate via <a href="#">LJM Library</a> Pico Technology PT-104 Platinum Resistance Data Logger, which communicate via <a href="#">PicoSDK/libusbpt104</a> TiePie (HS5, HS6, WS5) oscilloscopes, generators and I2C hosts, which communicate via <a href="#">LibTiePie SDK</a>

### 1.1.3 Sensor and Unit Conversion Utility

The Conversion Utility is a submodule that allows on the one hand a unified implementation of hardware-sensors and on the other hand provides a unified way to convert units. Furthermore it is possible to map two ranges on to each other. This can be useful to convert between for example and 4 - 20 mA and 0 - 10 V, both of them are common as sensor out- or input. Moreover, a subclass allows the mapping of a bit-range to any other range. For example a 12 bit number (0-4095) to 0 - 10. All utilities can be used with single numbers (`int`, `float`) as well as array-like structures containing single numbers (`np.array()`, `list`, `dict`, `tuple`).

Currently the following sensors are implemented:

- LEM LT 4000S
- LMT 70A

The following unit conversion classes are implemented:

- Temperature (Kelvin, Celsius, Fahrenheit)
- Pressure (Pascal, Bar, Atmosphere, Psi, Torr, Millimeter Mercury)

## 1.2 Documentation

Note: if you're planning to contribute to the `hvl_ccb` project read the **Contributing** section in the HVL CCB documentation.

Do either:

- read [HVL CCB documentation at RTD](#),

or

- build and read HVL CCB documentation locally; install first [Graphviz](#) (make sure to have the `dot` command in the executable search path) and the Python build requirements for documentation:

```
$ pip install docs/requirements.txt
```

and then either on Windows in Git BASH run:

```
$ ./make.sh docs
```

or from any other shell with GNU Make installed run:

```
$ make docs
```

The target index HTML ("`docs/_build/html/index.html`") should open automatically in your Web browser.

## 1.3 Credits

This package was created with [Cookiecutter](#) and the [audreyr/cookiecutter-pypackage](#) project template.



## **INSTALLATION**

### **2.1 Stable release**

To install HVL Common Code Base, run this command in your terminal:

```
$ pip install hvl_ccb
```

To install HVL Common Code Base with optional Python libraries that require manual installations of additional system libraries, you need to specify on installation extra requirements corresponding to these controllers. For instance, to install Python requirements for LabJack and TiePie devices, run:

```
$ pip install "hvl_ccb[tiepie,labjack]"
```

See below for the info about additional system libraries and the corresponding extra requirements.

To install all extra requirements run:

```
$ pip install "hvl_ccb[all]"
```

This is the preferred method to install HVL Common Code Base, as it will always install the most recent stable release.

If you don't have `pip` installed, this [Python installation guide](#) can guide you through the process.

### **2.2 From sources**

The sources for HVL Common Code Base can be downloaded from the [GitLab](#) repo.

You can either clone the repository:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

Or download the [tarball](#):

```
$ curl -OL https://gitlab.com/ethz_hvl/hvl_ccb/-/archive/master/hvl_ccb.tar.gz
```

Once you have a copy of the source, you can install it with:

```
$ pip install .
```

## 2.3 Additional system libraries

If you have installed *hvl\_ccb* with any of the extra features corresponding to device controllers, you must additionally install respective system library; these are:

Extra feature	Additional system library
labjack	<a href="#">LJM Library</a>
picotech	<a href="#">PicoSDK</a> (Windows) / <a href="#">libusbpt104</a> (Ubuntu/Debian)
tiepie	<a href="#">LibTiePie SDK</a>

For more details on installation of the libraries see docstrings of the corresponding `hvl_ccb` modules.

---

**CHAPTER  
THREE**

---

**USAGE**

To use HVL Common Code Base in a project:

```
import hvl_ccb
```



## API DOCUMENTATION


### 4.1 hvl\_ccb

#### 4.1.1 Subpackages

[hvl\\_ccb.comm](#)

**Submodules**

[hvl\\_ccb.comm.base](#)

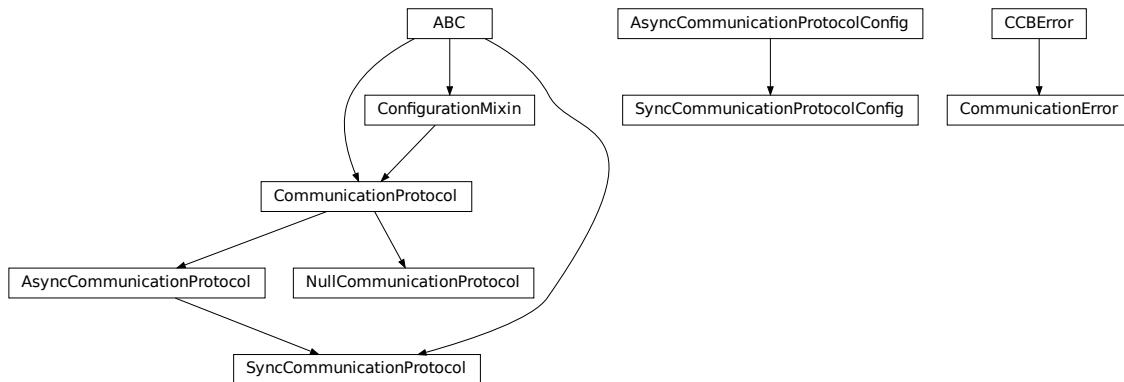

Module with base classes for communication protocols.

```
class AsyncCommunicationProtocol(config)
    Bases: CommunicationProtocol
    Abstract base class for asynchronous communication protocols
    static config_cls() → Type[AsyncCommunicationProtocolConfig]
        Return the default configdataclass class.

    Returns
        a reference to the default configdataclass class
```

**read()** → str

Read a single line of text as *str* from the communication.

**Returns**

text as *str* including the terminator, which can also be empty “”

**read\_all**(*n\_attempts\_max*: *Optional[int]* = *None*, *attempt\_interval\_sec*: *Optional[Union[int, float]]* = *None*) → *Optional[str]*

Read all lines of text from the connection till nothing is left to read.

**Parameters**

- **n\_attempts\_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt\_interval\_sec** – time between the reading attempts

**Returns**

A multi-line *str* including the terminator internally

**abstract read\_bytes()** → bytes

Read a single line as *bytes* from the communication.

This method uses *self.access\_lock* to ensure thread-safety.

**Returns**

a single line as *bytes* containing the terminator, which can also be empty b””

**read\_nonempty**(*n\_attempts\_max*: *Optional[int]* = *None*, *attempt\_interval\_sec*: *Optional[Union[int, float]]* = *None*) → *Optional[str]*

Try to read a non-empty single line of text as *str* from the communication. If the host does not reply or reply with white space only, it will return None.

**Returns**

a non-empty text as a *str* or *None* in case of an empty string

**Parameters**

- **n\_attempts\_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt\_interval\_sec** – time between the reading attempts

**read\_text()** → str

Read one line of text from the serial port. The input buffer may hold additional data afterwards, since only one line is read.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

**Returns**

String read from the serial port; “” if there was nothing to read.

**Raises**

[SerialCommunicationIOError](#) – when communication port is not opened

**read\_text\_nonempty**(*n\_attempts\_max*: *Optional[int]* = *None*, *attempt\_interval\_sec*: *Optional[Union[int, float]]* = *None*) → *Optional[str]*

Reads from the serial port, until a non-empty line is found, or the number of attempts is exceeded.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Attention: in contrast to *read\_text*, the returned answer will be stripped of a whitespace newline terminator at the end, if such terminator is set in the initial configuration (default).

**Parameters**

- **n\_attempts\_max** – maximum number of read attempts
- **attempt\_interval\_sec** – time between the reading attempts

**Returns**

String read from the serial port; ‘’ if number of attempts is exceeded or serial port is not opened.

**write(*text: str*)**

Write text as *str* to the communication.

**Parameters**

**text** – test as a *str* to be written

**abstract write\_bytes(*data: bytes*) → int**

Write data as *bytes* to the communication.

This method uses *self.access\_lock* to ensure thread-safety.

**Parameters**

**data** – data as *bytes*-string to be written

**Returns**

number of bytes written

**write\_text(*text: str*)**

Write text to the serial port. The text is encoded and terminated by the configured terminator.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

**Parameters**

**text** – Text to send to the port.

**Raises**

**SerialCommunicationIOError** – when communication port is not opened

**class AsyncCommunicationProtocolConfig(terminator: bytes = b'\n', encoding: str = 'utf-8', encoding\_error\_handling: str = 'strict', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10)**

Bases: **object**

Base configuration data class for asynchronous communication protocols

**clean\_values()****default\_n\_attempts\_read\_text\_nonempty: int = 10**

default number of attempts to read a non-empty text

**encoding: str = 'utf-8'**

Standard encoding of the connection. Typically this is **utf-8**, but can also be **latin-1** or something from here: <https://docs.python.org/3/library/codecs.html#standard-encodings>

**encoding\_error\_handling: str = 'strict'**

Encoding error handling scheme as defined here: <https://docs.python.org/3/library/codecs.html#error-handlers> By default strict error handling that raises **UnicodeError**.

**force\_value(*fieldname, value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

### Returns

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

### Returns

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

### Returns

a list of strings containing all required keys.

**terminator: bytes = b'\r\n'**

The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations.  
This defines the end of a single line.

**wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5**

time to wait between attempts of reading a non-empty text

## exception CommunicationError

Bases: [CCErrror](#)

**class CommunicationProtocol(config)**

Bases: [ConfigurationMixin](#), ABC

Communication protocol abstract base class.

Specifies the methods to implement for communication protocol, as well as implements some default settings and checks.

**access\_lock**

Access lock to use with context manager when accessing the communication protocol (thread safety)

**abstract close()**

Close the communication protocol

**abstract open()**

Open communication protocol

**class NullCommunicationProtocol(config)**

Bases: [CommunicationProtocol](#)

Communication protocol that does nothing.

**close()** → *None*

Void close function.

**static config\_cls()** → Type[*EmptyConfig*]

Empty configuration

**Returns**

*EmptyConfig*

**open()** → *None*

Void open function.

**class SyncCommunicationProtocol**(*config*)

Bases: *AsyncCommunicationProtocol*, ABC

Abstract base class for synchronous communication protocols with *query()*

**static config\_cls()** → Type[*SyncCommunicationProtocolConfig*]

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**query(command: str)** → Optional[str]

Send a command to the interface and handle the status message. Possibly raises an exception.

**Parameters**

**command** – Command to send


**Returns**

Answer from the interface, which can be None instead of an empty reply

**class SyncCommunicationProtocolConfig**(*terminator: bytes = b'\r\n'*, *encoding: str = 'utf-8'*,  
*encoding\_error\_handling: str = 'strict'*,  
*wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5*,  
*default\_n\_attempts\_read\_text\_nonempty: int = 10*)

Bases: *AsyncCommunicationProtocolConfig*

## hvl\_ccb.comm.labjack\_ljm

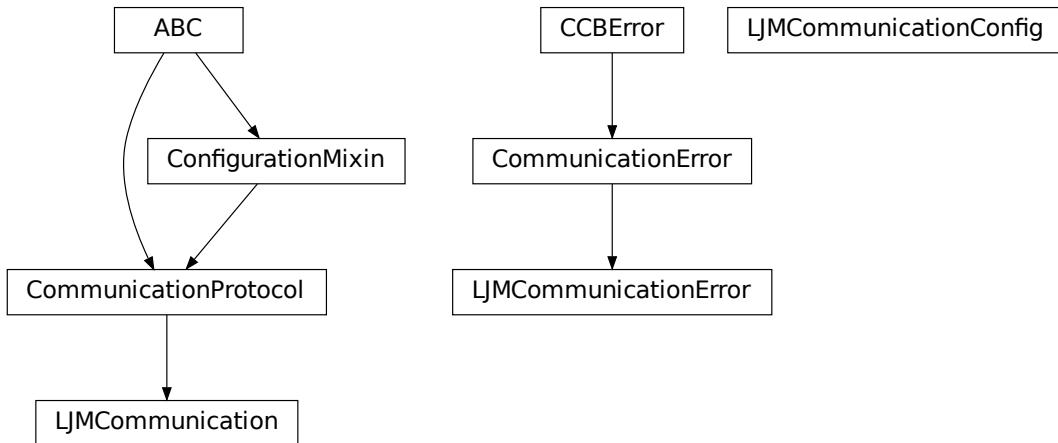

Communication protocol for LabJack using the LJM Library. Originally developed and tested for LabJack T7-PRO.

Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

**class LJMCommunication(configuration)**

Bases: `CommunicationProtocol`

Communication protocol implementing the LabJack LJM Library Python wrapper.

**close() → None**

Close the communication port.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**property is\_open: bool**

Flag indicating if the communication port is open.

**Returns**

*True* if the port is open, otherwise *False*

**open() → None**

Open the communication port.

**read\_name(\*names: str, return\_num\_type: ~typing.Type[~numbers.Real] = <class 'float'>) → Union[Real, Sequence[Real]]**

Read one or more input numeric values by name.

**Parameters**

- **names** – one or more names to read out from the LabJack

- **return\_num\_type** – optional numeric type specification for return values; by default *float*.

**Returns**

answer of the LabJack, either single number or multiple numbers in a sequence, respectively, when one or multiple names to read were given

**Raises**

**TypeError** – if read value of type not compatible with *return\_num\_type*

**write\_name**(*name*: str, *value*: Real) → *None*

Write one value to a named output.

**Parameters**

- **name** – String or with name of LabJack IO
- **value** – is the value to write to the named IO port

**write\_names**(*name\_value\_dict*: Dict[str, Real]) → *None*

Write more than one value at once to named outputs.

**Parameters**

**name\_value\_dict** – is a dictionary with string names of LabJack IO as keys and corresponding numeric values

**class LJMCommunicationConfig**(*device\_type*: Union[str, DeviceType] = 'ANY', *connection\_type*: Union[str, ConnectionType] = 'ANY', *identifier*: str = 'ANY')

Bases: *object*

Configuration dataclass for *LJMCommunication*.

**class ConnectionType**(*value*=<no\_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *AutoNumberNameEnum*

LabJack connection type.

**ANY** = 1

**ETHERNET** = 4

**TCP** = 3

**USB** = 2

**WIFI** = 5

**class DeviceType**(*value*=<no\_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *AutoNumberNameEnum*

LabJack device types.

Can be also looked up by ambiguous Product ID (*p\_id*) or by instance name: `python LabJackDeviceType(4)` is *LabJackDeviceType('T4')* `

**ANY** = 1

**T4** = 2

**T7** = 3

**T7\_PRO = 4**

**classmethod get\_by\_p\_id(p\_id: int) → Union[DeviceType, List[DeviceType]]**

Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

**Parameters**

**p\_id** – Product ID of a LabJack device

**Returns**

  Instance or list of instances of *LabJackDeviceType*

**Raises**

**ValueError** – when Product ID is unknown

**clean\_values() → None**

Performs value checks on device\_type and connection\_type.

**connection\_type: Union[str, ConnectionType] = 'ANY'**

Can be either string or of enum *ConnectionType*.

**device\_type: Union[str, DeviceType] = 'ANY'**

Can be either string ‘ANY’, ‘T7\_PRO’, ‘T7’, ‘T4’, or of enum *DeviceType*.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**identifier: str = 'ANY'**

The identifier specifies information for the connection to be used. This can be an IP address, serial number, or device name. See the LabJack docs (<https://labjack.com/support/software/api/ljm/function-reference/ljmopens/identifier-parameter>) for more information.

**is\_configdataclass = True**

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

  a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

  a list of strings containing all optional keys.

**classmethod required\_keys() → Sequence[str]**


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

  a list of strings containing all required keys.

**exception LJMCommunicationError**Bases: *CommunicationError*

Errors coming from LJMCommunication.

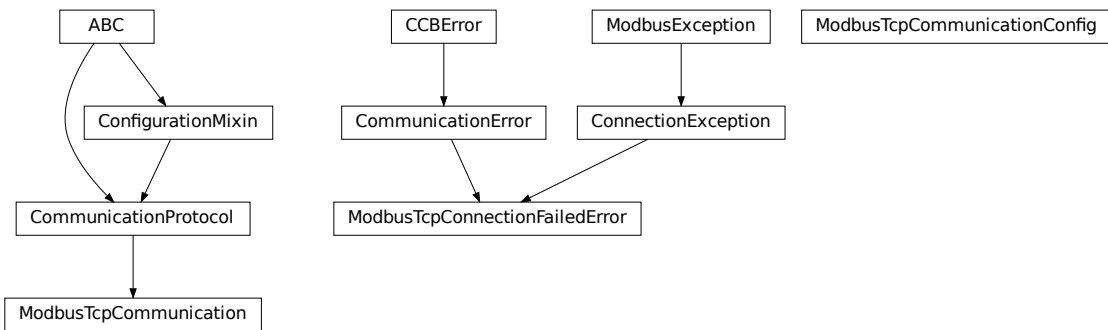**hvl\_ccb.comm.modbus\_tcp**Communication protocol for modbus TCP ports. Makes use of the `pymodbus` library.**class ModbusTcpCommunication(configuration)**Bases: *CommunicationProtocol*

Implements the Communication Protocol for modbus TCP.

**close()**

Close the Modbus TCP connection.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**open() → None**

Open the Modbus TCP connection.

**Raises**`ModbusTcpConnectionFailedError` – if the connection fails.**read\_holding\_registers(address: int, count: int) → List[int]**

Read specified number of register starting with given address and return the values from each register.

**Parameters**

- **address** – address of the first register
- **count** – count of registers to read

**Returns**list of `int` values**read\_input\_registers(address: int, count: int) → List[int]**

Read specified number of register starting with given address and return the values from each register in a list.

### Parameters

- **address** – address of the first register
- **count** – count of registers to read

### Returns

list of *int* values

**write\_registers**(*address*: *int*, *values*: *Union[List[int], int]*)

Write values from the specified address forward.

### Parameters

- **address** – address of the first register
- **values** – list with all values

**class ModbusTcpCommunicationConfig**(*host*: *Union[str, IPv4Address, IPv6Address]*, *unit*: *int*, *port*: *int* = 502)

Bases: *object*

Configuration dataclass for *ModbusTcpCommunication*.

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**host**: *Union[str, IPv4Address, IPv6Address]*

**is\_configdataclass** = *True*

**classmethod keys()** → *Sequence[str]*

Returns a list of all configdataclass fields key-names.

### Returns

a list of strings containing all keys.

**classmethod optional\_defaults()** → *Dict[str, object]*

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

### Returns

a list of strings containing all optional keys.

**port**: *int* = 502

**classmethod required\_keys()** → *Sequence[str]*


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

### Returns

a list of strings containing all required keys.

```
unit: int

exception ModbusTcpConnectionFailedError(string= "")
    Bases: ConnectionException, CommunicationError
    Error raised when the connection failed.
```

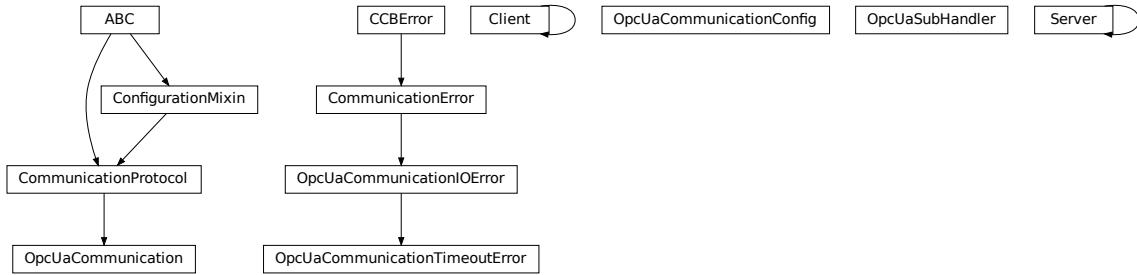**hvl\_ccb.comm.opc**

Communication protocol implementing an OPC UA connection. This protocol is used to interface with the “Cube” PLC from Siemens.

```
class Client(url: str, timeout: int = 4)
    Bases: Client
    disconnect()
    get_objects_node()
        Get Objects node of client. Returns a Node object.
    property is_open
    send_hello(*args, **kwargs)
class OpcUaCommunication(config)
    Bases: CommunicationProtocol
    Communication protocol implementing an OPC UA connection. Makes use of the package python-opcua.
    close() → None
        Close the connection to the OPC UA server.
    static config_cls()
        Return the default configdataclass class.

        Returns
            a reference to the default configdataclass class
    init_monitored_nodes(node_id: Union[object, Iterable], ns_index: int) → None
        Initialize monitored nodes.

        Parameters
            • node_id – one or more strings of node IDs; node IDs are always casted via str() method here, hence do not have to be strictly string objects.
```

- **ns\_index** – the namespace index the nodes belong to.

**Raises**

*OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

**property is\_open: bool**

Flag indicating if the communication port is open. —DEPRECATED! DO NOT USE!!!—

**Returns**

*True* if the port is open, otherwise *False*

**open() → None**

Open the communication to the OPC UA server.

**Raises**

*OpcUaCommunicationIOError* – when communication port cannot be opened.

**read(node\_id, ns\_index)**

Read a value from a node with id and namespace index.

**Parameters**

- **node\_id** – the ID of the node to read the value from
- **ns\_index** – the namespace index of the node

**Returns**

the value of the node object.

**Raises**

*OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

**write(node\_id, ns\_index, value) → None**

Write a value to a node with name **name**.

**Parameters**

- **node\_id** – the id of the node to write the value to.
- **ns\_index** – the namespace index of the node.
- **value** – the value to write.

**Raises**

*OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

```
class OpcUaCommunicationConfig(host: ~typing.Union[str, ~ipaddress.IPv4Address, ~ipaddress.IPv6Address],  
                               endpoint_name: str, port: int = 4840, sub_handler:  
                               ~hvl_ccb.comm.opc.OpcUaSubHandler =  
                               <hvl_ccb.comm.opc.OpcUaSubHandler object>, update_parameter:  
                               ~asyncua.ua.uaproto.Auto.CreateSubscriptionParameters =  
                               CreateSubscriptionParameters(RequestedPublishingInterval=1000,  
                               RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22,  
                               MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0),  
                               wait_timeout_retry_sec: ~typing.Union[int, float] = 1,  
                               max_timeout_retry_nr: int = 5)
```

Bases: **object**

Configuration dataclass for OPC UA Communciation.

**clean\_values()****endpoint\_name: str**

Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host: Union[str, IPv4Address, IPv6Address]**

Hostname or IP-Address of the OPC UA server.

**is\_configdataclass = True****classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**max\_timeout\_retry\_nr: int = 5**

Maximal number of call re-tries on underlying OPC UA client timeout error

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**port: int = 4840**

Port of the OPC UA server to connect to.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**sub\_handler: OpcUaSubHandler = <hvl\_ccb.comm.opc.OpcUaSubHandler object>**

object to use for handling subscriptions.

**update\_parameter: CreateSubscriptionParameters = CreateSubscriptionParameters(RequestedPublishingInterval=1000, RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22, MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0)**Values are given as a *ua.CreateSubscriptionParameters* as these parameters are requested by the OPC server. Other values will lead to an automatic revision of the parameters and a warning in the opc-logger, cf. MR !173

```

wait_timeout_retry_sec: Union[int, float] = 1
    Wait time between re-trying calls on underlying OPC UA client timeout error

exception OpcUaCommunicationIOError
    Bases: OSError, CommunicationError
    OPC-UA communication I/O error.

exception OpcUaCommunicationTimeoutError
    Bases: OpcUaCommunicationIOError
    OPC-UA communication timeout error.

class OpcUaSubHandler
    Bases: object
    Base class for subscription handling of OPC events and data change events. Override methods from this class to add own handling capabilities.


    To receive events from server for a subscription data_change and event methods are called directly from receiving thread. Do not do expensive, slow or network operation there. Create another thread if you need to do such a thing.

        datachange_notification(node, val, data)
        event_notification(event)

class Server(shelf_file=None, tloop=None)
    Bases: Server
    get_objects_node()
        Get Objects node of server. Returns a Node object.

```

## **hvl\_ccb.comm.serial**

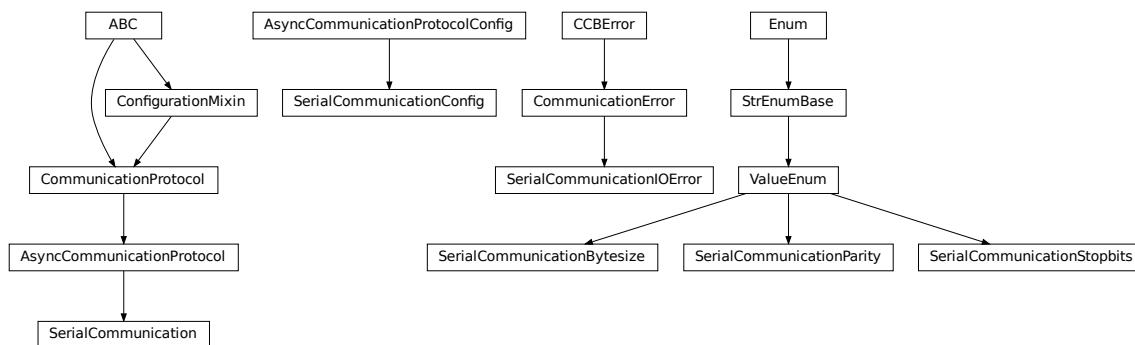

Communication protocol for serial ports. Makes use of the `pySerial` library.

```

class SerialCommunication(configuration)
    Bases: AsyncCommunicationProtocol
    Implements the Communication Protocol for serial ports.

```

**close()**

Close the serial connection.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**property is\_open: bool**

Flag indicating if the serial port is open.

**Returns**

*True* if the serial port is open, otherwise *False*

**open()**

Open the serial connection.

**Raises**

*SerialCommunicationIOError* – when communication port cannot be opened.

**read\_bytes() → bytes**

Read the bytes from the serial port till the terminator is found. The input buffer may hold additional lines afterwards.

This method uses *self.access\_lock* to ensure thread-safety.

**Returns**

Bytes read from the serial port; *b”* if there was nothing to read.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**read\_single\_bytes(size: int = 1) → bytes**

Read the specified number of bytes from the serial port. The input buffer may hold additional data afterwards.

**Returns**

Bytes read from the serial port; *b”* if there was nothing to read.

**write\_bytes(data: bytes) → int**

Write bytes to the serial port.

This method uses *self.access\_lock* to ensure thread-safety.

**Parameters**

**data** – data to write to the serial port

**Returns**

number of bytes written

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**class SerialCommunicationBytesize(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *ValueEnum*

Serial communication bytesize.

**EIGHTBITS = 8**

```
FIVEBITS = 5
SEVENBITS = 7
SIXBITS = 6

class SerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,
                                port: Optional[str] = None, baudrate: int = 9600, parity: Union[str,
                                SerialCommunicationParity] = SerialCommunicationParity.NONE,
                                stopbits: Union[int, float, SerialCommunicationStopbits] =
                                SerialCommunicationStopbits.ONE, bytesize: Union[int,
                                SerialCommunicationBytesize] =
                                SerialCommunicationBytesize.EIGHTBITS, timeout: Union[int, float] = 2)
```

Bases: [AsyncCommunicationProtocolConfig](#)

Configuration dataclass for [SerialCommunication](#).

#### **Bytesize**

alias of [SerialCommunicationBytesize](#)

#### **Parity**

alias of [SerialCommunicationParity](#)

#### **Stopbits**

alias of [SerialCommunicationStopbits](#)

#### **baudrate: int = 9600**

Baudrate of the serial port

#### **bytesize: Union[int, SerialCommunicationBytesize] = 8**

Size of a byte, 5 to 8

#### **clean\_values()**

#### **create\_serial\_port() → Serial**

Create a serial port instance according to specification in this configuration

#### **Returns**

Closed serial port instance

#### **force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### **Parameters**

- **fieldname** – name of the field
- **value** – value to assign

#### **classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

#### **Returns**

a list of strings containing all keys.

---

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

Parity to be used for the connection.

**port: Optional[str] = None**

Port is a string referring to a COM-port (e.g. 'COM3') or a URL. The full list of capabilities is found [on the pyserial documentation](#).

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**stopbits: Union[int, float, SerialCommunicationStopbits] = 1**

Stopbits setting, can be 1, 1.5 or 2.

**terminator\_str()** → str

**timeout: Union[int, float] = 2**

Timeout in seconds for the serial port

**exception SerialCommunicationIOError**

Bases: OSError, [CommunicationError](#)

Serial communication related I/O errors.

**class SerialCommunicationParity(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: [ValueEnum](#)

Serial communication parity.

**EVEN = 'E'**

**MARK = 'M'**

**NAMES = {'E': 'Even', 'M': 'Mark', 'N': 'None', 'O': 'Odd', 'S': 'Space'}**

**NONE = 'N'**

**ODD = 'O'**

**SPACE = 'S'**

**class SerialCommunicationStopbits(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**


Bases: [ValueEnum](#)

Serial communication stopbits.

**ONE = 1**

```
ONE_POINT_FIVE = 1.5  
TWO = 2
```

## hvl\_ccb.comm.tcp

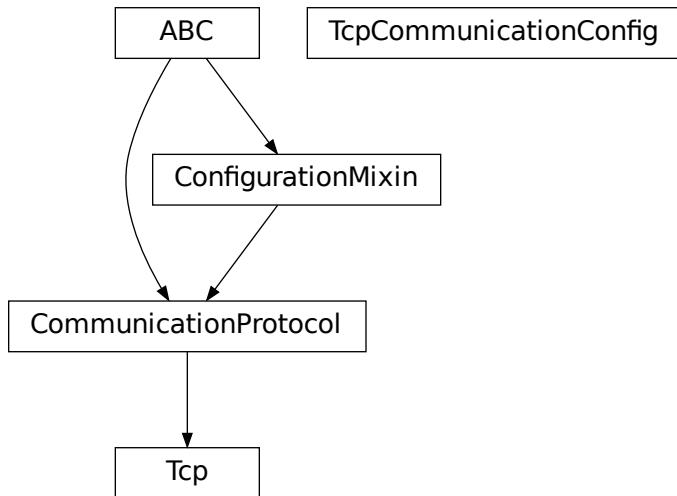

TCP communication protocol.

Makes use of the socket library.

```
class Tcp(configuration)  
    Bases: CommunicationProtocol  
    Tcp Communication Protocol.  
    close() → None  
        Close TCP connection.  
    static config_cls() → Type[TcpCommunicationConfig]  
        Return the default configdataclass class.  
        Returns  
            a reference to the default configdataclass class  
    open() → None  
        Open TCP connection.  
    read() → str  
        TCP read function :return: information read from TCP buffer formatted as string  
    write(command: str = "") → None  
        TCP write function :param command: command string to be sent :return: none  
class TcpCommunicationConfig(host: Union[str, IPv4Address, IPv6Address], port: int = 54321, bufsize: int = 1024)
```

Bases: `object`

Configuration dataclass for `TcpCommunication`.

**bufsize: int = 1024**

**clean\_values()**

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host: Union[str, IPv4Address, IPv6Address]**

**is\_configdataclass = True**

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**port: int = 54321**


**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

## hvl\_ccb.comm.telnet

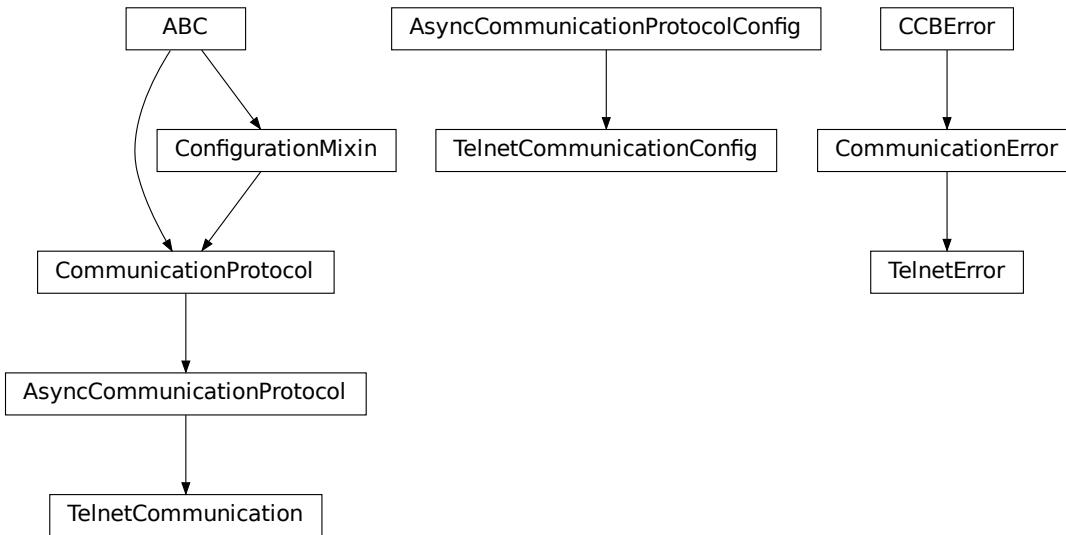

Communication protocol for telnet. Makes use of the `telnetlib` library.

**class TelnetCommunication(configuration)**

Bases: `AsyncCommunicationProtocol`

Implements the Communication Protocol for telnet.

**close()**

Close the telnet connection unless it is not closed.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**property is\_open: bool**

Is the connection open?

**Returns**

True for an open connection

**open()**

Open the telnet connection unless it is not yet opened.

**read\_bytes() → bytes**

Read data as `bytes` from the telnet connection.

**Returns**

data from telnet connection

**Raises**

`TelnetError` – when connection is not open, raises an Error during the communication

**write\_bytes**(*data: bytes*)

Write the data as *bytes* to the telnet connection.

**Parameters**

- **data** – Data to be sent.

**Raises**

- **TelnetError** – when connection is not open, raises an Error during the communication

```
class TelnetCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                    encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                    Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,
                                    host: Optional[Union[str, IPv4Address, IPv6Address]] = None, port: int =
                                    0, timeout: Union[int, float] = 0.2)
```

Bases: *AsyncCommunicationProtocolConfig*

Configuration dataclass for *TelnetCommunication*.

**clean\_values()****create\_telnet()** → *Optional[Telnet]*

Create a telnet client :return: Opened Telnet object or None if connection is not possible

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host: Optional[Union[str, IPv4Address, IPv6Address]] = None**

Host to connect to can be localhost or

**classmethod keys()** → *Sequence[str]*

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → *Dict[str, object]*

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**port: int = 0**

Port at which the host is listening

**classmethod required\_keys()** → *Sequence[str]*

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.


```
timeout: Union[int, float] = 0.2
```

Timeout for reading a line

**exception TelnetError**

Bases: OSError, *CommunicationError*

Telnet communication related errors.

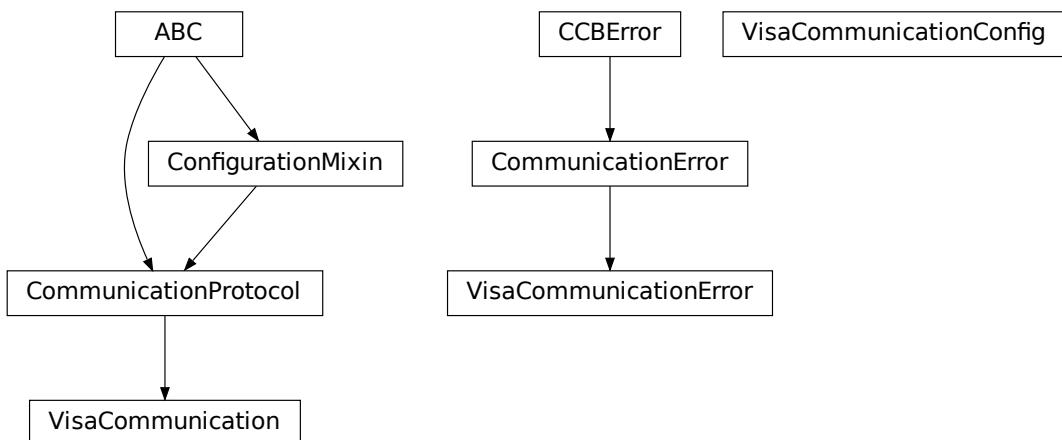**hvl\_ccb.comm.visa**

Communication protocol for VISA. Makes use of the pyvisa library. The backend can be NI-Visa or pyvisa-py.

Information on how to install a VISA backend can be found here: [https://pyvisa.readthedocs.io/en/master/getting\\_nivisa.html](https://pyvisa.readthedocs.io/en/master/getting_nivisa.html)

So far only TCPIP SOCKET and TCPIP INSTR interfaces are supported.

**class VisaCommunication(configuration)**

Bases: *CommunicationProtocol*

Implements the Communication Protocol for VISA / SCPI.

**MULTI\_COMMANDS\_MAX = 5**

The maximum of commands that can be sent in one round is 5 according to the VISA standard.

**MULTI\_COMMANDS\_SEPARATOR = ';'**

The character to separate two commands is ; according to the VISA standard.

**WAIT\_AFTER\_WRITE = 0.08**

Small pause in seconds to wait after write operations, allowing devices to really do what we tell them before continuing with further tasks.

**close() → None**

Close the VISA connection and invalidates the handle.

**static config\_cls()** → Type[*VisaCommunicationConfig*]

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**open()** → *None*

Open the VISA connection and create the resource.

**query(\*commands: str)** → Union[str, Tuple[str, ...]]

A combination of write(message) and read.

**Parameters**

**commands** – list of commands

**Returns**

list of values

**Raises**

**VisaCommunicationError** – when connection was not started, or when trying to issue too many commands at once.

**spoll()** → int

Execute serial poll on the device. Reads the status byte register STB. This is a fast function that can be executed periodically in a polling fashion.

**Returns**

integer representation of the status byte

**Raises**

**VisaCommunicationError** – when connection was not started

**write(\*commands: str)** → *None*

Write commands. No answer is read or expected.

**Parameters**

**commands** – one or more commands to send

**Raises**

**VisaCommunicationError** – when connection was not started

**class VisaCommunicationConfig(host: Union[str, IPv4Address, IPv6Address], interface\_type: Union[str, InterfaceType], board: int = 0, port: int = 5025, timeout: int = 5000, chunk\_size: int = 204800, open\_timeout: int = 1000, write\_termination: str = '\n', read\_termination: str = '\n', visa\_backend: str = '')**

Bases: object

*VisaCommunication* configuration dataclass.

**class InterfaceType(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *AutoNumberNameEnum*

Supported VISA Interface types.

**TCPIP\_INSTR = 2**

VXI-11 protocol

**TCPIP\_SOCKET = 1**

VISA-Raw protocol

**address**(*host: str, port: Optional[int] = None, board: Optional[int] = None*) → str

Address string specific to the VISA interface type.

**Parameters**

- **host** – host IP address
- **port** – optional TCP port
- **board** – optional board number

**Returns**

address string

**property address: str**

Address string depending on the VISA protocol's configuration.

**Returns**

address string corresponding to current configuration

**board: int = 0**

Board number is typically 0 and comes from old bus systems.

**chunk\_size: int = 204800**

Chunk size is the allocated memory for read operations. The standard is 20kB, and is increased per default here to 200kB. It is specified in bytes.

**clean\_values()**

**force\_value**(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host: Union[str, IPv4Address, IPv6Address]**

**interface\_type: Union[str, InterfaceType]**

Interface type of the VISA connection, being one of *InterfaceType*.

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**open\_timeout: int = 1000**

Timeout for opening the connection, in milli seconds.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

```
port: int = 5025
    TCP port, standard is 5025.

read_termination: str = '\n'
    Read termination character.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns
    a list of strings containing all required keys.

timeout: int = 5000
    Timeout for commands in milli seconds.

visa_backend: str = ''
    Specifies the path to the library to be used with PyVISA as a backend. Defaults to None, which is NI-VISA (if installed), or pyvisa-py (if NI-VISA is not found). To force the use of pyvisa-py, specify '@py' here.

write_termination: str = '\n'
    Write termination character.

exception VisaCommunicationError
    Bases: OSError, CommunicationError
    Base class for VisaCommunication errors.
```

## Module contents

Communication protocols subpackage.


### **hvl\_ccb.dev**

#### **Subpackages**

##### **hvl\_ccb.dev.cube**

#### **Submodules**

## `hvl_ccb.dev(cube).base`

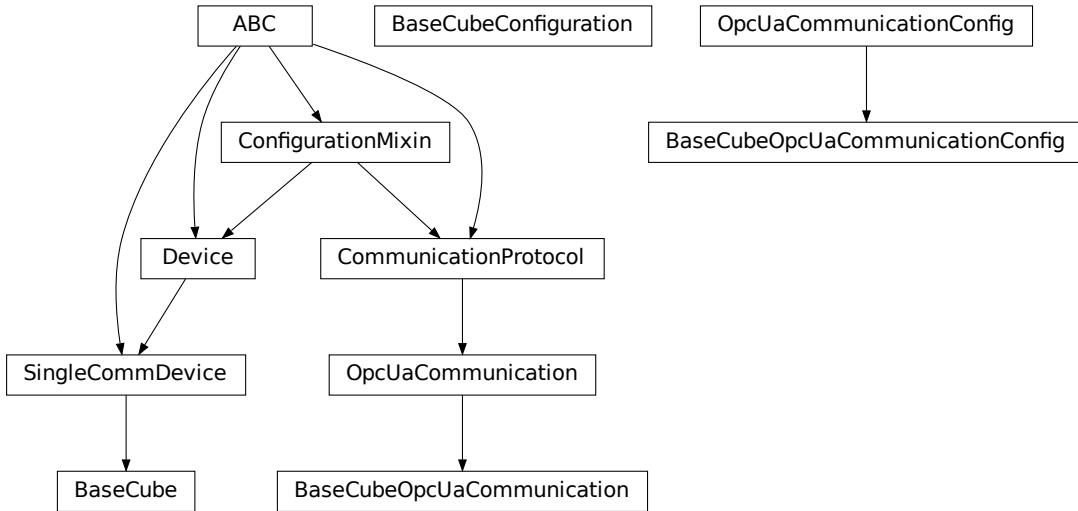

Classes for the BaseCube device.

`class BaseCube(com, dev_config=None)`

Bases: `SingleCommDevice`

Base class for Cube variants.

`OPC_MAX_YEAR = 2089`

`OPC_MIN_YEAR = 1990`

`active_alarms(human_readable: bool = True) → List[Union[int, str]]`

Displays all active alarms / messages.

**Parameters**

`human_readable` – `True` for human readable message, `False` for corresponding integer

**Returns**

list with active alarms

`property breakdown_detection_active: bool`

Get the state of the breakdown detection functionality. Returns True if it is enabled, False otherwise.

**Returns**

state of the breakdown detection functionality

`breakdown_detection_reset() → None`

Reset the breakdown detection circuitry so that it is ready to detect breakdowns again.

`property breakdown_detection_triggered: bool`

See if breakdown detection unit has been triggered. Returns True if it is triggered, False otherwise.

**Returns**

trigger status of the breakdown detection unit

**property cee16\_socket**

Read the on-state of the IEC CEE16 three-phase power socket.

**Returns**

the on-state of the CEE16 power socket

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**classmethod datetime\_to\_opc(*time\_dt*: datetime) → List[int]**

Converts python datetime format into opc format (list of 8 integers) as defined in the following link: <https://support.industry.siemens.com/cs/mdm/109798671?c=133950752267&lc=de-WW> Each byte corresponds to one list entry. [yy, MM, dd, hh, mm, ss, milliseconds, weekday] Milliseconds and Weekday are not used, as this precision / information is not needed. The conversion of the numbers is special. Each decimal number is treated as it would be a hex-number and then converted back to decimal. This is tested with the used PLC in the BaseCube. yy: 0 to 99 (0 -> 2000, 89 -> 2089, 90 -> 1990, 99 -> 1999) MM: 1 to 12 dd: 1 to 31 hh: 0 to 23 mm: 0 to 59 ss: 0 to 59

**Parameters**

**time\_dt** – time to be converted

**Returns**

time in opc list format

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**display\_message\_board() → None**

Display 15 newest messages

**display\_status\_board() → None**

Display status board.

**door\_1\_status**

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

**door\_2\_status**

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

**door\_3\_status**

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

**earthing\_rod\_1\_status**

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

**earthing\_rod\_2\_status**

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

**earthing\_rod\_3\_status**

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

**property operate: Optional[bool]**

Indicates if ‘operate’ is activated. ‘operate’ means locket safety circuit, red lamps, high voltage on and locked safety switches.

**Returns**

*True* if operate is activated (RED\_OPERATE), *False* if ready is deactivated (RED\_READY),  
*None* otherwise

**quit\_error()** → *None*

Quits errors that are active on the Cube.

**read(node\_id: str)**

Local wrapper for the OPC UA communication protocol read method.

**Parameters**

**node\_id** – the id of the node to read.

**Returns**

the value of the variable

**property ready: Optional[bool]**

Indicates if ‘ready’ is activated. ‘ready’ means locket safety circuit, red lamps, but high voltage still off.

**Returns**

*True* if ready is activated (RED\_READY), *False* if ready is deactivated (GREEN\_READY),  
*None* otherwise

**set\_message\_board(msgs: List[str], display\_board: bool = True)** → *None*

Fills messages into message board that display that 15 newest messages with a timestamp.

**Parameters**

- **msgs** – list of strings
- **display\_board** – display 15 newest messages if *True* (default)

**Raises**

**ValueError** – if there are too many messages or the positions indices are invalid.

**set\_status\_board(msgs: List[str], pos: Optional[List[int]] = None, clear\_board: bool = True, display\_board: bool = True)** → *None*

Sets and displays a status board. The messages and the position of the message can be defined.

**Parameters**

- **msgs** – list of strings
- **pos** – list of integers [0...14]
- **clear\_board** – clear unspecified lines if *True* (default), keep otherwise
- **display\_board** – display new status board if *True* (default)

**Raises**

**ValueError** – if there are too many messages or the positions indices are invalid.

**start()** → *None*

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

**property status: SafetyStatus**

Get the safety circuit status of the Cube. This methods is for the user.

**Returns**

the safety status of the Cube's state machine.

**stop() → None**

Stop the Cube device. Deactivates the remote control and closes the communication protocol.

**Raises**

**CubeStopError** – when the cube is not in the correct status to stop the operation

**t13\_socket\_1**

Set and get the state of a SEV T13 power socket.

**t13\_socket\_2**

Set and get the state of a SEV T13 power socket.

**t13\_socket\_3**

Set and get the state of a SEV T13 power socket.

**write(node\_id, value) → None**

Local wrapper for the OPC UA communication protocol write method.

**Parameters**

- **node\_id** – the id of the node to write
- **value** – the value to write to the variable

```
class BaseCubeConfiguration(namespace_index: int = 3, polling_delay_sec: Union[int, float] = 5.0,
                            polling_interval_sec: Union[int, float] = 1.0, timeout_status_change:
                            Union[int, float] = 6, timeout_interval: Union[int, float] = 0.1,
                            noise_level_measurement_channel_1: Union[int, float] = 100,
                            noise_level_measurement_channel_2: Union[int, float] = 100,
                            noise_level_measurement_channel_3: Union[int, float] = 100,
                            noise_level_measurement_channel_4: Union[int, float] = 100)
```

Bases: object

Configuration dataclass for the BaseCube devices.

**clean\_values()****force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True****classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

```
namespace_index: int = 3
```

Namespace of the OPC variables, typically this is 3 (coming from Siemens)

```
noise_level_measurement_channel_1: Union[int, float] = 100
```

```
noise_level_measurement_channel_2: Union[int, float] = 100
```

```
noise_level_measurement_channel_3: Union[int, float] = 100
```

```
noise_level_measurement_channel_4: Union[int, float] = 100
```

```
classmethod optional_defaults() → Dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

```
polling_delay_sec: Union[int, float] = 5.0
```

```
polling_interval_sec: Union[int, float] = 1.0
```

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

```
timeout_interval: Union[int, float] = 0.1
```

```
timeout_status_change: Union[int, float] = 6
```

```
class BaseCubeOpcUaCommunication(config)
```

Bases: *OpcUaCommunication*

Communication protocol specification for BaseCube devices.

```
static config_cls()
```

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
class BaseCubeOpcUaCommunicationConfig(host: ~typing.Union[str, ~ipaddress.IPv4Address,
~ipaddress.IPv6Address], endpoint_name:
~hvl_ccb.dev(cube.constants._CubeOpcEndpoint =
_CubeOpcEndpoint.BASE_CUBE, port: int = 4840, sub_handler:
~hvl_ccb.comm.opc.OpcUaSubHandler =
<hvl_ccb.dev(cube.base._BaseCubeSubscriptionHandler object>,
update_parameter:
~asyncua.ua.uaproto.protocol_auto.CreateSubscriptionParameters =
CreateSubscriptionParameters(RequestedPublishingInterval=1000,
RequestedLifetimeCount=300,
RequestedMaxKeepAliveCount=22,
MaxNotificationsPerPublish=10000, PublishingEnabled=True,
Priority=0), wait_timeout_retry_sec: ~typing.Union[int, float] =
1, max_timeout_retry_nr: int = 5)
```

Bases: *OpcUaCommunicationConfig*

Communication protocol configuration for OPC UA, specifications for the BaseCube devices.

**endpoint\_name:** *\_CubeOpcEndpoint* = 'BaseCube'

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

#### Returns

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### Returns

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


#### Returns

a list of strings containing all required keys.

**sub\_handler:** *OpcUaSubHandler* = <hvl\_ccb.dev(cube.base.\_BaseCubeSubscriptionHandler object>

Subscription handler for data change events

## hvl\_ccb.dev(cube.constants

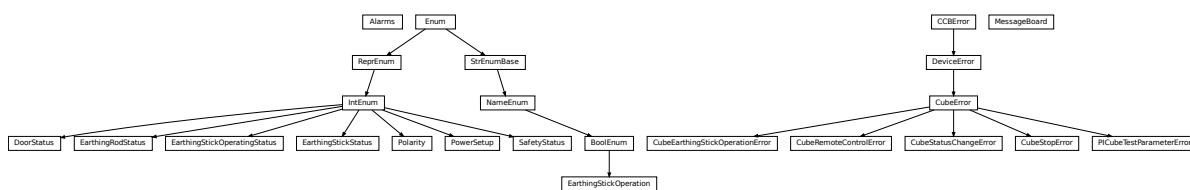

Constants, variable names for the BaseCube OPC-connected devices.

**exception CubeEarthlingStickOperationError**

Bases: *CubeError*

**exception CubeError**

Bases: *DeviceError*

**exception CubeRemoteControlError**

Bases: *CubeError*

**exception CubeStatusChangeError**

Bases: *CubeError*

**exception CubeStopError**

Bases: *CubeError*

**class DoorStatus**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *IntEnum*

Possible status values for doors.

**CLOSED = 2**

Door is closed, but not locked.

**ERROR = 4**

Door has an error or was opened in locked state (either with emergency stop or from the inside).

**INACTIVE = 0**

not enabled in BaseCube HMI setup, this door is not supervised.

**LOCKED = 3**

Door is closed and locked (safe state).

**OPEN = 1**

Door is open.

**class EarthingRodStatus**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *IntEnum*

Possible status values for earthing rods.

**EXPERIMENT\_BLOCKED = 0**

earthing rod is somewhere in the experiment and blocks the start of the experiment

**EXPERIMENT\_READY = 1**

earthing rod is hanging next to the door, experiment is ready to operate

**class EarthingStickOperatingStatus**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *IntEnum*

Operating Status for an earthing stick. Stick can be used in auto or manual mode.

**AUTO = 0**

**MANUAL = 1**

**class EarthingStickOperation**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *BoolEnum*

Operation of the earthing stick in manual operating mode. Can be closed or opened.

```

CLOSE = True
OPEN = False

class EarthingStickStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                           start=1, boundary=None)
Bases: IntEnum
Status of an earthing stick. These are the possible values in the status integer e.g. in _EarthingStick.status.

CLOSED = 1
ERROR = 3
INACTIVE = 0
OPEN = 2

class MessageBoard(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                     boundary=None)
Bases: _LineEnumBase
Variable NodeID strings for message board lines.

LINE_1 = '"DB_OPC_Connection"."Is_status_Line_1"'
LINE_10 = '"DB_OPC_Connection"."Is_status_Line_10"'
LINE_11 = '"DB_OPC_Connection"."Is_status_Line_11"'
LINE_12 = '"DB_OPC_Connection"."Is_status_Line_12"'
LINE_13 = '"DB_OPC_Connection"."Is_status_Line_13"'
LINE_14 = '"DB_OPC_Connection"."Is_status_Line_14"'
LINE_15 = '"DB_OPC_Connection"."Is_status_Line_15"'
LINE_2 = '"DB_OPC_Connection"."Is_status_Line_2"'
LINE_3 = '"DB_OPC_Connection"."Is_status_Line_3"'
LINE_4 = '"DB_OPC_Connection"."Is_status_Line_4"'
LINE_5 = '"DB_OPC_Connection"."Is_status_Line_5"'
LINE_6 = '"DB_OPC_Connection"."Is_status_Line_6"'
LINE_7 = '"DB_OPC_Connection"."Is_status_Line_7"'
LINE_8 = '"DB_OPC_Connection"."Is_status_Line_8"'
LINE_9 = '"DB_OPC_Connection"."Is_status_Line_9"'

exception PICubeTestParameterError
Bases: CubeError

```

```
class Polarity(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
               boundary=None)
```

Bases: IntEnum

An enumeration.

**NEGATIVE = 0**

**POSITIVE = 1**

```
class PowerSetup(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
```

Bases: IntEnum

Possible power setups corresponding to the value of variable Power.setup. The values for slope\_min are experimentally defined, below these values the slope is more like a staircase

The name of the first argument needs to be ‘value’, otherwise the IntEnum is not working correctly.

**AC\_100KV = 3**

**AC\_150KV = 4**

**AC\_200KV = 5**

**AC\_50KV = 2**

**DC\_140KV = 7**

**DC\_280KV = 8**

**EXTERNAL\_SOURCE = 1**

**IMPULSE\_140KV = 9**

**NO\_SOURCE = 0**

**POWER\_INVERTER\_220V = 6**

```
STOP_SAFETY_STATUSES: Tuple[SafetyStatus, ...] = (<SafetyStatus.GREEN\_NOT\_READY: 1>,
                                                       <SafetyStatus.GREEN\_READY: 2>)
```

BaseCube’s safety statuses required to close the connection to the device.

```
class SafetyStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                     boundary=None)
```

Bases: IntEnum

Safety status values that are possible states returned from [hvl\\_ccb.dev.cube.base.BaseCube.status\(\)](#). These values correspond to the states of the BaseCube’s safety circuit statemachine.

**ERROR = 6**

**GREEN\_NOT\_READY = 1**

**GREEN\_READY = 2**


**INITIALIZING = 0**

**QUICK\_STOP = 5**

```
RED_OPERATE = 4
```

```
RED_READY = 3
```

## hvl\_ccb.dev(cube.picube

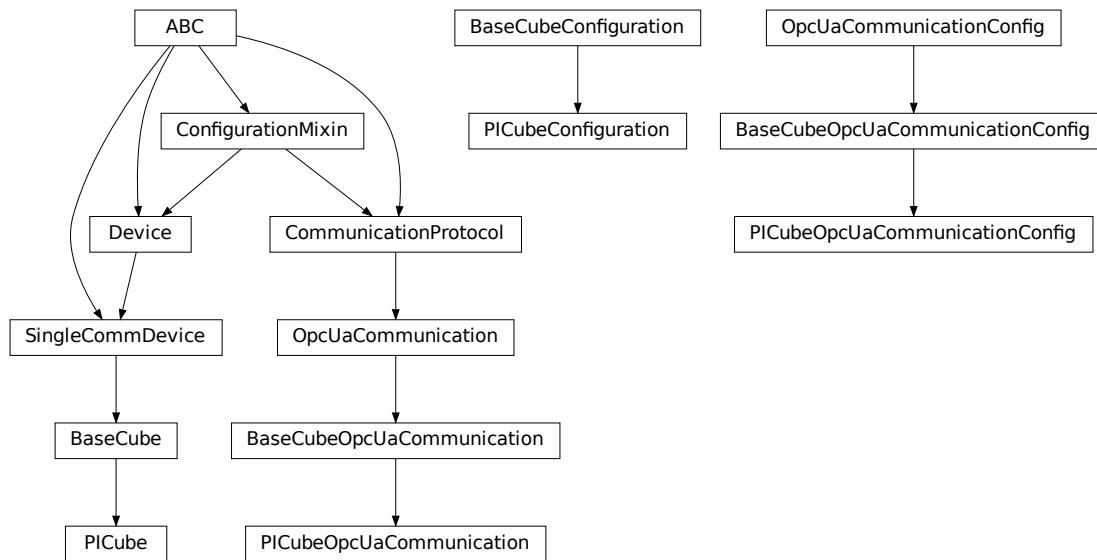

A PICube is a BaseCube with build in Power Inverter

```
class PICube(com, dev_config=None)
```

Bases: *BaseCube*

Variant of the BaseCube with build in Power Inverter

```
static config_cls()
```

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
property current_primary: float
```

Read the current primary current at the output of the frequency converter (before transformer).

**Returns**

primary current in A

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

```
property frequency: float
```

Read the electrical frequency of the current PICube setup.

**Returns**

the frequency in Hz

**property operate: Optional[bool]**

Indicates if ‘operate’ is activated. ‘operate’ means locket safety circuit, red lamps, high voltage on and locked safety switches.

**Returns**

*True* if operate is activated (RED\_OPERATE), *False* if ready is deactivated (RED\_READY),  
*None* otherwise

**property polarity: Optional[Polarity]**

Polarity of a DC setup. :return: if a DC setup is programmed the polarity is returned, else None.

**property power\_setup: PowerSetup**

Return the power setup selected in the PICube’s settings.

**Returns**

the power setup

**property voltage\_actual: float**

Reads the actual measured voltage and returns the value in V.

**Returns**

the actual voltage of the setup in V.

**property voltage\_max: float**

Reads the maximum voltage of the setup and returns in V.

**Returns**

the maximum voltage of the setup in V.

**property voltage\_primary: float**

Read the current primary voltage at the output of the frequency converter (before transformer).

**Returns**

primary voltage in V

```
class PICubeConfiguration(namespace_index: int = 3, polling_delay_sec: Union[int, float] = 5.0,
                           polling_interval_sec: Union[int, float] = 1.0, timeout_status_change: Union[int, float] = 6, timeout_interval: Union[int, float] = 0.1,
                           noise_level_measurement_channel_1: Union[int, float] = 100, noise_level_measurement_channel_2: Union[int, float] = 100,
                           noise_level_measurement_channel_3: Union[int, float] = 100, noise_level_measurement_channel_4: Union[int, float] = 100,
                           timeout_test_parameters: 'Number' = 2.0)
```

Bases: *BaseCubeConfiguration*

**clean\_values()****force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**timeout\_test\_parameters:** Union[int, float] = 2.0

**class PICubeOpcUaCommunication(config)**

Bases: *BaseCubeOpcUaCommunication*

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class PICubeOpcUaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], endpoint\_name: '\_CubeOpcEndpoint' = <\_CubeOpcEndpoint.PI\_CUBE: 'PICube'>, port: int = 4840, sub\_handler: hvl\_ccb.comm.opc.OpcUaSubHandler = <hvl\_ccb.dev.cube.base.\_BaseCubeSubscriptionHandler object at 0x7f883ea90f90>, update\_parameter: asynccua.ua.uaproto.protocol\_auto.CreateSubscriptionParameters = CreateSubscriptionParameters(RequestedPublishingInterval=1000, RequestedLifetimeCount=300, RequestedMaxKeepAliveCount=22, MaxNotificationsPerPublish=10000, PublishingEnabled=True, Priority=0), wait\_timeout\_retry\_sec: Union[int, float] = 1, max\_timeout\_retry\_nr: int = 5)**

Bases: *BaseCubeOpcUaCommunicationConfig*

**endpoint\_name: \_CubeOpcEndpoint = 'PICube'**

Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.


## Module contents

Cube package with implementation for system versions from 2019 on (new concept with hard-PLC Siemens S7-1500 as CPU).

## hvl\_ccb.dev.fluke884x

### Submodules

#### hvl\_ccb.dev.fluke884x.base

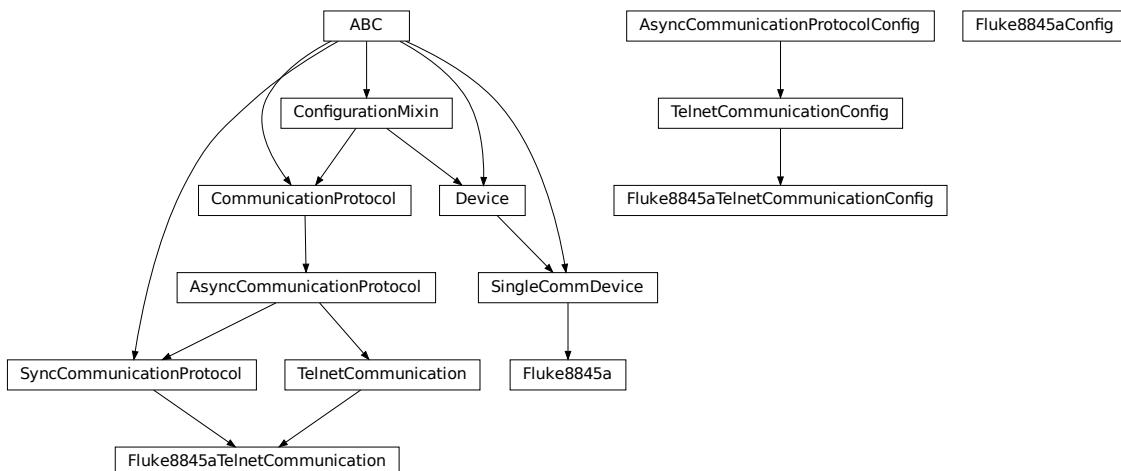

Python module for the Fluke8845a Multimeter. The communication to the device is through Telnet. 8845A/8846A Programmers Manual is available in the following link. All page numbers mentioned in this script refer to this manual. [https://download.flukecal.com/pub/literature/8845A\\_pmeng0300.pdf](https://download.flukecal.com/pub/literature/8845A_pmeng0300.pdf)

```
class Fluke8845a(com, dev_config=None)
```

Bases: *SingleCommDevice*

Device class to control Fluke8845a

```
DISPLAY_MAX_LENGTH = 12
```

```
ac_current_range
```

```
ac_voltage_range
```

```
activate_remote_mode() → None
```

Page 66

Places the Meter in the remote mode for RS-232 or Ethernet remote control. All front-panel keys, except the local key, are disabled.

```
clear_display_message() → None
```

Page 59

Clears the displayed message on the Meter's display.

```
clear_error_queue() → None
```

Page 62

Sets all bits to zero in the Meter's status byte register and all event registers. Also clears the error queue

```
static config_cls() → Type[Fluke8845aConfig]
```

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
current_filter
```

```
dc_current_range
```

```
dc_voltage_range
```

```
static default_com_cls() → Type[Fluke8845aTelnetCommunication]
```

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

```
property display_enable: bool
```

Page 59

get if the display is enabled or not fluke answer string “1” for ON and “0” for off bool(int(“1”)) = 1 and  
bool(int(“0”)) = 0

**Returns**

bool enabled = True, else False

```
property display_message: str
```

Page 59

Retrieves the text sent to the Meter's display.

**fetch()** → float

Page 36

Transfer stored readings to output buffer

**four\_wire\_resistance\_range**

**frequency\_aperture**

**property identification: str**

Page 60

Queries “\*IDN?” and returns the identification string of the connected device.

**Returns**

the identification string of the connected device e.g. “FLUKE, 8845A, 2540017, 08/02/10-11:53”

**initiate\_trigger()** → *None*

Set trigger system to wait-for-trigger

**measure()** → float

Page 42

Taking measurement

Once the Meter has been configured for a measurement, the INITiate command causes the Meter to take a measurement when the trigger condition have been met. To process readings from the Meter’s internal memory to the output buffer, send the Meter a FETCh? command.

**property measurement\_function: MeasurementFunction**

input\_function getter, query what the input function is

**Raises**

*Fluke8845aUnknownCommandError* – if the input function is unknown

**period\_aperture**

**reset()** → *None*

Page 60

resets the meter to its power-up configuration

**start()** → *None*

Start this device as recommended by the manual

**stop()** → *None*

Stop this device. Disables access and closes the communication protocol.

**trigger()** → *None*

Causes the meter to trigger a measurement when paused

**property trigger\_delay: int**

input\_trigger\_delay getter, query what the input trigger delay is in second answer format from Fluke: string, ‘+1.0000000E+00’, so convert to float and then to int

**Returns**

input trigger delay in second

```

property trigger_source: TriggerSource
    input_trigger_source getter, query what the input trigger source is

Raises
    Fluke8845aUnknownCommandError – if the input trigger source is unknown

two_wire_resistance_range

voltage_filter

class Fluke8845aConfig(name: str = 'Fluke 1')
    Bases: object
    Config for Fluke8845a
    name: the name of the device
    clean_values()
        Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.
    force_value(fieldname, value)
        Forces a value to a dataclass field despite the class being frozen.
        NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.
    Parameters
        • fieldname – name of the field
        • value – value to assign
    is_configdataclass = True
    classmethod keys() → Sequence[str]
        Returns a list of all configdataclass fields key-names.

    Returns
        a list of strings containing all keys.

    name: str = 'Fluke 1'

    classmethod optional_defaults() → Dict[str, object]
        Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns
        a list of strings containing all optional keys.

    classmethod required_keys() → Sequence[str]
        Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns
        a list of strings containing all required keys.

class Fluke8845aTelnetCommunication(configuration)
    Bases: TelnetCommunication, SyncCommunicationProtocol

```

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**query(command: str) → str**

Send a command to the interface and handle the status message. Eventually raises an error.

**Parameters**

**command** – Command to send

**Raises**

**Fluke8845aError** – if the connection is broken

**Returns**

Answer from the interface

**class Fluke8845aTelnetCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8', encoding\_error\_handling: str = 'strict', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10, host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address, NoneType] = None, port: int = 3490, timeout: Union[int, float] = 0.2)**

Bases: *TelnetCommunicationConfig*

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**port: int = 3490**

Port at which Fluke 8845a is listening

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


**Returns**

a list of strings containing all required keys.

```
terminator: bytes = b'\r'
```

The terminator is CR

## `hvl_ccb.dev.fluke884x.constants`

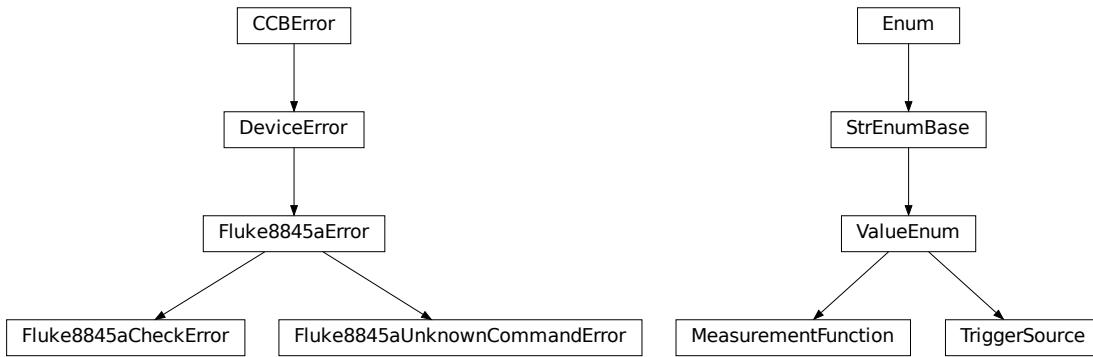

Constants, ValueEnum: MeasurementFunction and TriggerSoruce Descriptors for range, filter and aperture

### `exception Fluke8845aCheckError`

Bases: `Fluke8845aError`

### `exception Fluke8845aError`

Bases: `DeviceError`

### `exception Fluke8845aUnknownCommandError`

Bases: `Fluke8845aError`

```
class MeasurementFunction(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                           start=1, boundary=None)
```

Bases: `ValueEnum`

Page 40

Sets the Meter function. This command must be followed by the INIT and FETCh? commands to cause the meter to take a measurement.

`CURRENT_AC = 'CURR:AC'`

`CURRENT_DC = 'CURR'`

`DIODE = 'DIOD'`

`FOUR_WIRE_RESISTANCE = 'FRES'`

`FREQUENCY = 'FREQ'`

`PERIOD = 'PER'`

`TWO_WIRE_RESISTANCE = 'RES'`

`VOLTAGE_AC = 'VOLT:AC'`

```
VOLTAGE_DC = 'VOLT'
```

```
class TriggerSource(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                     boundary=None)
```

Bases: [ValueEnum](#)

Page 57

**BUS:** Sets the Meter to expect a trigger through the IEEE-488 bus or upon execution of a \*TRG command

**IMM:** Selects Meter's internal triggering system


**EXT:** Sets the Meter to sense triggers through the trigger jack on the rear panel of the Meter

```
BUS = 'BUS'
```

```
EXTERNAL = 'EXT'
```

```
IMMEDIATE = 'IMM'
```

## [hvl\\_ccb.dev.fluke884x.ranges](#)

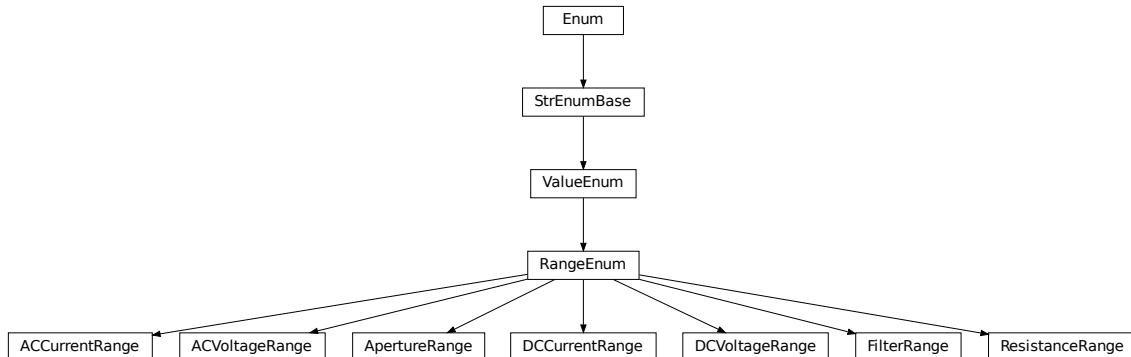

Ranges, RangeEnum for Fluke8845a device

```
class ACCurrentRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                      boundary=None)
```

Bases: [RangeEnum](#)

possible measurement ranges for AC current with unit Ampere

```
FOUR_HUNDRED_MILLI_AMPERE = 0.4
```

```
ONE_AMPERE = 1.0
```

```
ONE_HUNDRED_MILLI_AMPERE = 0.1
```

```
TEN_AMPERE = 10.0
```

```
TEN_MILLI_AMPERE = 0.01
```

```
THREE_AMPERE = 3.0
```

```
class ACVoltageRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
boundary=None)
```

Bases: [RangeEnum](#)

possible measurement ranges for AC voltage with unit volt

```
HUNDRED_VOLT = 100.0
```

```
ONE_HUNDRED_MILLI_VOLT = 0.1
```

```
ONE_VOLT = 1.0
```

```
SEVEN_HUNDRED_FIFTY_VOLT = 750.0
```

```
TEN_VOLT = 10.0
```

```
class ApertureRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
boundary=None)
```

Bases: [RangeEnum](#)

Page 46

Sets the gate time for the frequency/period function to the value  
 10ms = 4 1/2 digits  
 100ms = 5 1/2 digits  
 1s = 6 1/2 digits

```
ONE_HUNDRED_MILLI_SECOND = 0.1
```

```
ONE_SECOND = 1.0
```

```
TEN_MILLI_SECOND = 0.01
```

```
class DCCurrentRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
boundary=None)
```

Bases: [RangeEnum](#)

possible measurement ranges for DC current with unit Ampere

```
FOUR_HUNDRED_MILLI_AMPERE = 0.4
```

```
ONE_AMPERE = 1.0
```

```
ONE_HUNDRED_MICRO_AMPERE = 0.0001
```

```
ONE_HUNDRED_MILLI_AMPERE = 0.1
```

```
ONE_MILLI_AMPERE = 0.001
```

```
TEN_AMPERE = 10.0
```

```
TEN_MILLI_AMPERE = 0.01
```

```
THREE_AMPERE = 3.0
```

```
class DCVoltageRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                     boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for DC voltage with unit volt

**HUNDRED\_VOLT = 100.0**

**ONE\_HUNDRED\_MILLI\_VOLT = 0.1**

**ONE\_THOUSAND\_VOLT = 1000.0**

**ONE\_VOLT = 1.0**

**TEN\_VOLT = 10.0**

```
class FilterRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
```

Bases: *RangeEnum*

Page 47

```
Sets the appropriate filter for the frequency specified by <n>
High pass filter
For `VOLTAGE_AC`: <n> Hz to 300 kHz
For `CURRENT_AC`: <n> Hz to 10 kHz
parameters <n> = 3 slow filter
                  20 medium filter
                  200 fast filter
For `CURRENT_AC` and `VOLTAGE_AC`
```

**FAST\_FILTER = 200.0**

**MEDIUM\_FILTER = 20.0**

**SLOW\_FILTER = 3.0**

```
class ResistanceRange(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                      boundary=None)
```

Bases: *RangeEnum*

possible measurement ranges for resistance with unit Ohm

**ONE\_HUNDRED\_MILLION\_OHM = 1000000000.0**

**ONE\_HUNDRED\_OHM = 100.0**

**ONE\_HUNDRED\_THOUSAND\_OHM = 100000.0**

**ONE\_MILLION\_OHM = 1000000.0**

**ONE\_THOUSAND\_OHM = 1000.0**

**TEN\_MILLION\_OHM = 10000000.0**

**TEN\_THOUSAND\_OHM = 10000.0**


## Module contents

Fluke 8845A multimeter implementation using Telnet communication

### [hvl\\_ccb.dev.highland\\_t560](#)

#### Submodules

##### [hvl\\_ccb.dev.highland\\_t560.base](#)

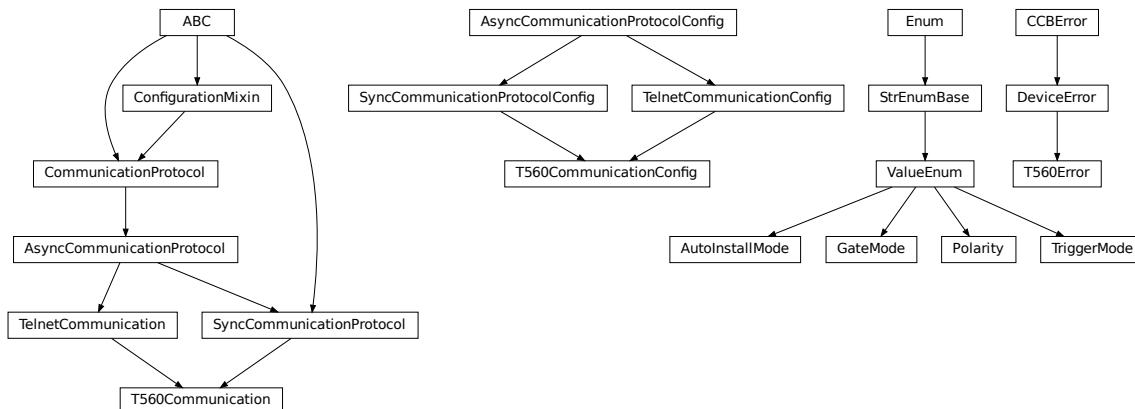

Module containing base device and communication classes and enums.

Communication with device is performed via its ethernet port and a Telnet connection.

```
class AutoInstallMode(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                      boundary=None)
```

Bases: [ValueEnum](#)

Modes for installing configuration settings to the device.

**INSTALL** = 1

**OFF** = 0

**QUEUE** = 2

```
class GateMode(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
               boundary=None)
```

Bases: [ValueEnum](#)

Available T560 gate modes

**INPUT** = 'INP'

**OFF** = 'OFF'

**OUTPUT** = 'OUT'

```
class Polarity(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
               boundary=None)
```

Bases: [ValueEnum](#)

Possible channel polarity states

```
ACTIVE_HIGH = 'POS'
```

```
ACTIVE_LOW = 'NEG'
```

```
class T560Communication(configuration)
```

Bases: [SyncCommunicationProtocol](#), [TelnetCommunication](#)

Communication class for T560. It uses a TelnetCommunication with the SyncCommunicationProtocol

```
static config_cls()
```

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
query(command: str) → str
```

Send a command to the device and handle the response.

For device setting queries, response will be ‘OK’ if successful, or ‘?’ if setting cannot be carried out, raising an error.

**Parameters**

**command** – Command string to be sent

**Raises**

[T560Error](#) – if no response is received, or if the device responds with an error message.

**Returns**

Response from the device.

```
class T560CommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8', encoding_error_handling:
                                str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                default_n_attempts_read_text_nonempty: int = 10, host: Union[str,
                                ipaddress.IPv4Address, ipaddress.IPv6Address, NoneType] = None, port: int
                                = 2000, timeout: Union[int, float] = 0.2)
```

Bases: [SyncCommunicationProtocolConfig](#), [TelnetCommunicationConfig](#)

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**port: int = 2000**

Port at which the host is listening

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**terminator: bytes = b'\r'**

The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations.  
This defines the end of a single line.

**exception T560Error**

Bases: *DeviceError*

T560 related errors.

**class TriggerMode(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *ValueEnum*

Available T560 trigger modes

**COMMAND = 'REM'**

**EXT\_FALLING\_EDGE = 'NEG'**


**EXT\_RISING\_EDGE = 'POS'**

**INT\_SYNTHESIZER = 'SYN'**

**OFF = 'OFF'**

**hvl\_ccb.dev.highland\_t560.channel**

Module for controlling pulse output channels A, B, C and D.

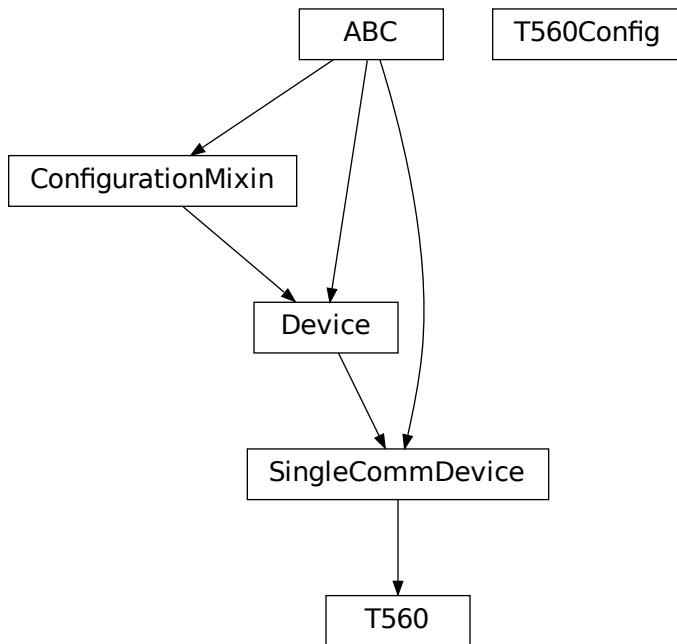**hvl\_ccb.dev.highland\_t560.device**

Module for controlling device, including TRIG, CLOCK and GATE I/Os.

```

class T560(com, dev_config=None)
    Bases: SingleCommDevice

    activate_clock_output()
        Outputs 10 MHz clock signal

    property auto_install_mode: AutoInstallMode
        Check the autoinstall settings of the T560. The autoinstall mode sets how changes to device settings are applied. See manual section 4.7.2 for more information about these modes.

    property ch_a: _Channel
        Channel A of T560

    property ch_b: _Channel
        Channel B of T560

    property ch_c: _Channel
        Channel C of T560

    property ch_d: _Channel
        Channel D of T560

    static config_cls()
        Return the default configdataclass class.
  
```

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**disarm\_trigger()**

Disarm DDG by disabling all trigger sources.

**fire\_trigger()**

Fire a software trigger.

**property frequency: float**

The frequency of the timing cycle in Hz.

**property gate\_mode: GateMode**

Check the mode setting of the GATE I/O port.

**property gate\_polarity: Polarity**

Check the polarity setting of the GATE I/O port.

**load\_device\_configuration()**

Load the settings saved in nonvolatile memory.

**property period: float**

The period of the timing cycle (time between triggers) in seconds.

**save\_device\_configuration()**

Save the current settings to nonvolatile memory.

**property trigger\_level**

Get external trigger level.

**property trigger\_mode**

Get device trigger source.

**use\_external\_clock()**

Finds and accepts an external clock signal to the CLOCK input

**class T560Config**

Bases: object

**auto\_install\_mode = 1****clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**is\_configdataclass** = True

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.


## Module contents

This module establishes methods for interfacing with the Highland Technology T560-2 via its ethernet adapter with a telnet communication protocol.

The T560 is a small digital delay & pulse generator. It outputs up to four individually timed pulses with 10-ps precision, given an internal or external trigger.

This module introduces methods for configuring channels, gating, and triggering. Further documentation and a more extensive command list may be obtained from:

<https://www.highlandtechnology.com/DSS/T560DS.shtml>

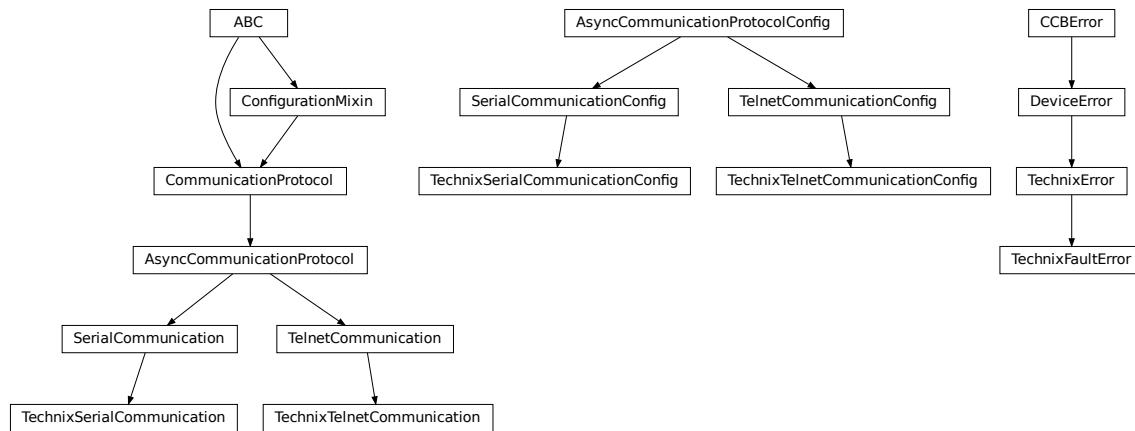**hvl\_ccb.dev.technix****Submodules****hvl\_ccb.dev.technix.base**

Communication and auxiliary classes for Technix

**exception TechnixError**

Bases: `DeviceError`

Technix related errors.

**exception TechnixFaultError**

Bases: `TechnixError`

Raised when the fault flag was detected while the interlock is closed

**class TechnixSerialCommunication(configuration)**

Bases: `_TechnixCommunication`, `SerialCommunication`

Serial communication for Technix

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
class TechnixSerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, port:
                                         Optional[str] = None, baudrate: int = 9600, parity: Union[str,
                                         SerialCommunicationParity] =
                                         SerialCommunicationParity.NONE, stopbits: Union[int, float,
                                         SerialCommunicationStopbits] =
                                         SerialCommunicationStopbits.ONE, bytesize: Union[int,
                                         SerialCommunicationBytesize] =
                                         SerialCommunicationBytesize.EIGHTBITS, timeout: Union[int,
                                         float] = 2)
```

Bases: `_TechnixCommunicationConfig`, `SerialCommunicationConfig`

Configuration for the serial communication for Technix

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**class TechnixTelnetCommunication**(*configuration*)

Bases: `TelnetCommunication`, `_TechnixCommunication`

Telnet communication for Technix

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
class TechnixTelnetCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, host:
                                         Optional[Union[str, IPv4Address, IPv6Address]] = None, port:
                                         int = 4660, timeout: Union[int, float] = 0.2)
```

Bases: `_TechnixCommunicationConfig`, `TelnetCommunicationConfig`

Configuration for the telnet communication for Technix

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

#### Returns

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### Returns

a list of strings containing all optional keys.

**port: int = 4660**


Port at which Technix is listening

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

#### Returns

a list of strings containing all required keys.

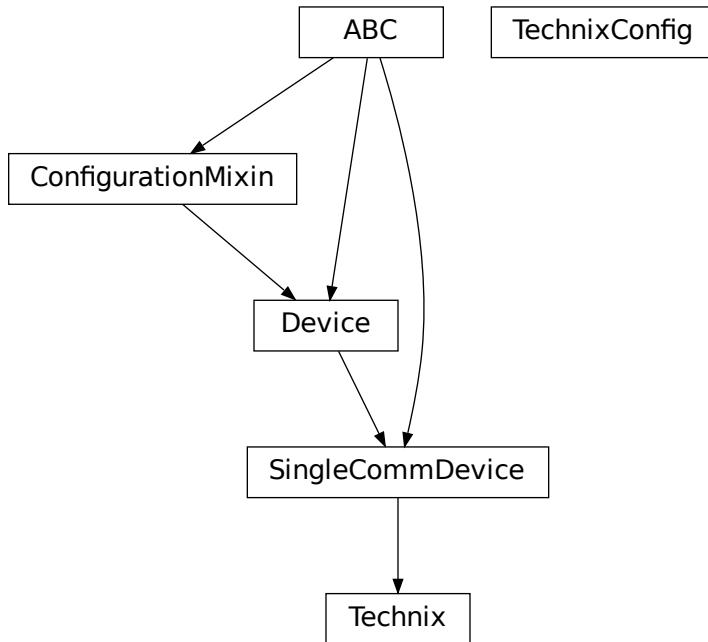**hvl\_ccb.dev.technix.device**

its corresponding configuration class

**class Technix(com, dev\_config)**

Bases: *SingleCommDevice*

Device class to control capacitor chargers from Technix

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**property current: Optional[Union[int, float]]**

Actual current of the output in A

**default\_com\_cls() → Union[Type[TechnixSerialCommunication], Type[TechnixTelnetCommunication]]**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**property inhibit: Optional[bool]**

Is the output of the voltage inhibited? The output stage can still be active.

**property is\_started: bool**

Is the device started?

The device class *Technix* and

```
property max_current: Union[int, float]
    Maximal output current of the hardware in A
property max_voltage: Union[int, float]
    Maximal output voltage of the hardware in V
property open_interlock: Optional[bool]
    Is the interlock open? (in safe mode)
property output: Optional[bool]
    State of the high voltage output
query_status(*, _retry: bool = False)
    Query the status of the device.

Returns
    This function returns nothing

property remote: Optional[bool]
    Is the device in remote control mode?

start()
    Start the device and set it into the remote controllable mode. The high voltage is turn off, and the status poller is started.

property status: Optional[_Status]
    The status of the device with the different states as sub-fields

stop()
    Stop the device. The status poller is stopped and the high voltage output is turn off.

property voltage: Optional[Union[int, float]]
    Actual voltage at the output in V
property voltage_regulation: Optional[bool]
    Status if the output is in voltage regulation mode (or current regulation)

class TechnixConfig(communication_channel:
    Union[Type[hvl_ccb.dev.technix.base.TechnixSerialCommunication],
          Type[hvl_ccb.dev.technix.base.TechnixTelnetCommunication]], max_voltage: Union[int, float], max_current: Union[int, float], polling_interval_sec: Union[int, float] = 4, post_stop_pause_sec: Union[int, float] = 1, register_pulse_time: Union[int, float] = 0.1, read_output_while_polling: bool = False)

Bases: object

clean_values()
    Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

communication_channel: Union[Type[TechnixSerialCommunication], Type[TechnixTelnetCommunication]]
    communication channel between computer and Technix

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.
```

### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass** = True

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

### Returns

a list of strings containing all keys.

**max\_current: Union[int, float]**

Maximal Output current

**max\_voltage: Union[int, float]**

Maximal Output voltage

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

### Returns

a list of strings containing all optional keys.

**polling\_interval\_sec: Union[int, float] = 4**

Polling interval in s to maintain to watchdog of the device

**post\_stop\_pause\_sec: Union[int, float] = 1**

Time to wait after stopping the device

**read\_output\_while\_polling: bool = False**

Read output voltage and current within the polling event

**register\_pulse\_time: Union[int, float] = 0.1**

Time for pulsing a register

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

### Returns

a list of strings containing all required keys.

## Module contents

Device classes for “RS 232” and “Ethernet” interfaces, which are used to control power supplies from Technix. Manufacturer homepage: <https://www.technix-hv.com>

The regulated power supplies series and capacitor chargers series from Technix are series of low and high voltage direct current power supplies as well as capacitor chargers. The class *Technix* is tested with a CCR10KV-7,5KJ via an ethernet connection as well as a CCR15-P-2500-OP via a serial connection. Check the code carefully before using it with other devices or device series

This Python package may support the following interfaces from Technix:

- Remote Interface RS232

- Ethernet Remote Interface
- Optic Fiber Remote Interface

## **hvl\_ccb.dev.tiepie**

### **Submodules**

#### **hvl\_ccb.dev.tiepie.base**

**NOTE:** `libtiepie` tries to load the `libtiepie` SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match `docs/hvl_ccb.dev.tiepie.XXX.rst` and remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.base
    :parts: 1

.. automodule:: hvl_ccb.dev.tiepie.base
    :members:
    :undoc-members:
    :show-inheritance:
```

#### **hvl\_ccb.dev.tiepie.channel**

**NOTE:** `libtiepie` tries to load the `libtiepie` SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match `docs/hvl_ccb.dev.tiepie.XXX.rst` and remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.channel
    :parts: 1

.. automodule:: hvl_ccb.dev.tiepie.channel
    :members:
    :undoc-members:
    :show-inheritance:
```

### **hvl\_ccb.dev.tiepie.device**

**NOTE:** `libtiepie` tries to load the `libtiepie` SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match `docs/hvl_ccb.dev.tiepie.XXX.rst` and remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.device
:parts: 1

.. automodule:: hvl_ccb.dev.tiepie.device
:members:
:undoc-members:
:show-inheritance:
```

### **hvl\_ccb.dev.tiepie.generator**

**NOTE:** `libtiepie` tries to load the `libtiepie` SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match `docs/hvl_ccb.dev.tiepie.XXX.rst` and remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.generator
:parts: 1

.. automodule:: hvl_ccb.dev.tiepie.generator
:members:
:undoc-members:
:show-inheritance:
```

### **hvl\_ccb.dev.tiepie.i2c**

**NOTE:** `libtiepie` tries to load the `libtiepie` SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match `docs/hvl_ccb.dev.tiepie.XXX.rst` and remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.i2c
:parts: 1

.. automodule:: hvl_ccb.dev.tiepie.i2c
:members:
:undoc-members:
:show-inheritance:
```

## hvl\_ccb.dev.tiepie.oscilloscope

**NOTE:** libtiepie tries to load the libtiepie SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match docs/hvl\_ccb.dev.tiepie.XXX.rst and remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.oscilloscope
:parts: 1

.. automodule:: hvl_ccb.dev.tiepie.oscilloscope
:members:
:undoc-members:
:show-inheritance:
```

## hvl\_ccb.dev.tiepie.utils

**NOTE:** libtiepie tries to load the libtiepie SDK during import. Thus, the API docs can only be generated in a system with the latter installed.

To build the API docs for this subpackage locally edit all files that match docs/hvl\_ccb.dev.tiepie.XXX.rst and remove the `.. code-block::` directive preceding the following directives:


```
.. inheritance-diagram:: hvl_ccb.dev.tiepie.utils
:parts: 1

.. automodule:: hvl_ccb.dev.tiepie.utils
:members:
:undoc-members:
:show-inheritance:
```

## Module contents

### Submodules

## `hvl_ccb.dev.base`

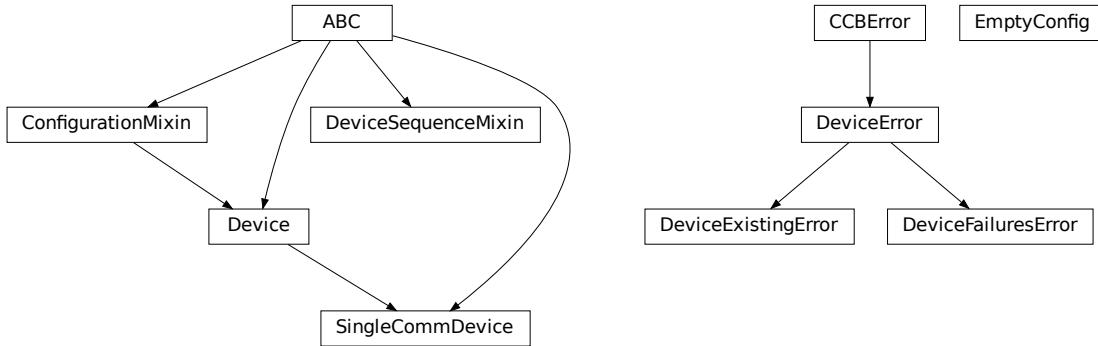

Module with base classes for devices.

**class Device(*dev\_config=None*)**

Bases: *ConfigurationMixin*, ABC

Base class for devices. Implement this class for a concrete device, such as measurement equipment or voltage sources.

Specifies the methods to implement for a device.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**abstract start() → *None***

Start or restart this Device. To be implemented in the subclass.

**abstract stop() → *None***

Stop this Device. To be implemented in the subclass.

**exception DeviceError**

Bases: *CCBError*

**exception DeviceExistingError**

Bases: *DeviceError*

Error to indicate that a device with that name already exists.

**exception DeviceFailuresError(*failures: Dict[str, Exception]*, \*args)**

Bases: *DeviceError*

Error to indicate that one or several devices failed.

**failures: Dict[str, Exception]**

A dictionary of named devices failures (exceptions).

**class DeviceSequenceMixin(*devices*: Dict[str, Device])**

Bases: ABC

Mixin that can be used on a device or other classes to provide facilities for handling multiple devices in a sequence.

**add\_device(*name*: str, *device*: Device) → None**

Add a new device to the device sequence.

**Parameters**

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

**Raises**

*DeviceExistingError* –

**devices\_failed\_start: Dict[str, Device]**

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices\_failed\_stop*.

**devices\_failed\_stop: Dict[str, Device]**

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices\_failed\_start*.

**get\_device(*name*: str) → Device**

Get a device by name.

**Parameters**

- **name** – is the name of the device.

**Returns**

the device object from this sequence.

**get\_devices() → List[Tuple[str, Device]]**

Get list of name, device pairs according to current sequence.

**Returns**

A list of tuples with name and device each.

**remove\_device(*name*: str) → Device**

Remove a device from this sequence and return the device object.

**Parameters**

- **name** – is the name of the device.

**Returns**

device object or *None* if such device was not in the sequence.

**Raises**

**ValueError** – when device with given name was not found

**start() → None**

Start all devices in this sequence in their added order.

**Raises**

*DeviceFailuresError* – if one or several devices failed to start

**stop()** → *None*

Stop all devices in this sequence in their reverse order.

**Raises**

*DeviceFailuresError* – if one or several devices failed to stop

**class EmptyConfig**

Bases: `object`

Empty configuration dataclass that is the default configuration for a Device.

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**class SingleCommDevice(*com*, *dev\_config=None*)**

Bases: `Device`, ABC

Base class for devices with a single communication protocol.

**property com**

Get the communication protocol of this device.

**Returns**

an instance of CommunicationProtocol subtype

**abstract static default\_com\_cls()** → Type[*CommunicationProtocol*]

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device


**start()** → *None*

Open the associated communication protocol.

**stop()** → *None*

Close the associated communication protocol.

## hvl\_ccb.dev.crylas

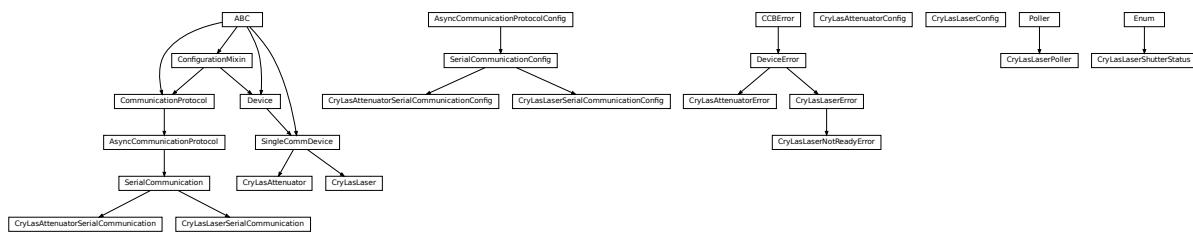

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: [https://www.crylas.de/products/pulsed\\_laser.html](https://www.crylas.de/products/pulsed_laser.html)

**class CryLasAttenuator(*com, dev\_config=None*)**

Bases: *SingleCommDevice*

Device class for the CryLas laser attenuator.

**property attenuation: Union[int, float]**

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**set\_attenuation(*percent: Union[int, float]*)** → *None*

Set the percentage of attenuated light (inverse of `set_transmission`). :param percent: percentage of attenuation, number between 0 and 100 :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

**set\_init\_attenuation()**

Sets the attenuation to its configured initial/default value

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**set\_transmission(percent: Union[int, float]) → None**

Set the percentage of transmitted light (inverse of set\_attenuation). :param percent: percentage of transmitted light :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

**start() → None**

Open the com, apply the config value ‘init\_attenuation’

**Raises**

*SerialCommunicationIOError* – when communication port cannot be opened

**property transmission: Union[int, float]**

**class CryLasAttenuatorConfig(init\_attenuation: Union[int, float] = 0, response\_sleep\_time: Union[int, float] = 1)**

Bases: object

Device configuration dataclass for CryLas attenuator.

**clean\_values()**

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**init\_attenuation: Union[int, float] = 0**

**is\_configdataclass = True**

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**response\_sleep\_time:** Union[int, float] = 1

**exception CryLasAttenuatorError**

Bases: *DeviceError*

General error with the CryLas Attenuator.

**class CryLasAttenuatorSerialCommunication(configuration)**

Bases: *SerialCommunication*

Specific communication protocol implementation for the CryLas attenuator. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class CryLasAttenuatorSerialCommunicationConfig(terminator: bytes = b'', encoding: str = 'utf-8', encoding\_error\_handling: str = 'strict', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)**

Bases: *SerialCommunicationConfig*

**baudrate: int = 9600**

Baudrate for CryLas attenuator is 9600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

CryLas attenuator does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**stopbits: Union[int, SerialCommunicationStopbits] = 1**

CryLas attenuator uses one stop bit

**terminator: bytes = b''**

No terminator

**timeout: Union[int, float] = 3**

use 3 seconds timeout as default

**class CryLasLaser(com, dev\_config=None)**

Bases: *SingleCommDevice*

CryLas laser controller device class.

**class AnswersShutter(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *Enum*

Standard answers of the CryLas laser controller to ‘Shutter’ command passed via *com*.

**CLOSED = 'Shutter inaktiv'**

**OPENED = 'Shutter aktiv'**

**class AnswersStatus(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *Enum*

Standard answers of the CryLas laser controller to ‘STATUS’ command passed via *com*.

**ACTIVE = 'STATUS: Laser active'**

**HEAD = 'STATUS: Head ok'**

**INACTIVE = 'STATUS: Laser inactive'**

**READY = 'STATUS: System ready'**

**TEC1 = 'STATUS: TEC1 Regulation ok'**

```

TEC2 = 'STATUS: TEC2 Regulation ok'

class LaserStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
    Bases: Enum
    Status of the CryLas laser

    READY_ACTIVE = 2
    READY_INACTIVE = 1
    UNREADY_INACTIVE = 0

    property is_inactive
    property is_ready

class RepetitionRates(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                      start=1, boundary=None)
    Bases: IntEnum
    Repetition rates for the internal software trigger in Hz

    HARDWARE = 0
    SOFTWARE_INTERNAL_SIXTY = 60
    SOFTWARE_INTERNAL_TEN = 10
    SOFTWARE_INTERNAL_TWENTY = 20

ShutterStatus
    alias of CryLasLaserShutterStatus

close_shutter() → None
    Close the laser shutter.

Raises
    • SerialCommunicationIOError – when communication port is not opened
    • CryLasLaserError – if success is not confirmed by the device

static config_cls()
    Return the default configdataclass class.

    Returns
        a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns
        the type of the standard communication protocol for this device

get_pulse_energy_and_rate() → Tuple[int, int]
    Use the debug mode, return the measured pulse energy and rate.

    Returns
        (energy in micro joule, rate in Hz)

```

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not answer the query

**laser\_off()** → *None*

Turn the laser off.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**laser\_on()** → *None*

Turn the laser on.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserNotReadyError* – if the laser is not ready to be turned on
- *CryLasLaserError* – if success is not confirmed by the device

**open\_shutter()** → *None*

Open the laser shutter.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**set\_init\_shutter\_status()** → *None*

Open or close the shutter, to match the configured shutter\_status.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**set\_pulse\_energy(energy: int)** → *None*

Sets the energy of pulses (works only with external hardware trigger). Proceed with small energy steps, or the regulation may fail.

**Parameters**

**energy** – energy in micro joule

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not confirm success

**set\_repetition\_rate(rate: Union[int, RepetitionRates])** → *None*

Sets the repetition rate of the internal software trigger.

**Parameters**

**rate** – frequency (Hz) as an integer

**Raises**

- *ValueError* – if rate is not an accepted value in RepetitionRates Enum

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

**start()** → *None*

Opens the communication protocol and configures the device.

**Raises**

- **SerialCommunicationIOError** – when communication port cannot be opened

**stop()** → *None*

Stops the device and closes the communication protocol.

**Raises**

- **SerialCommunicationIOError** – if com port is closed unexpectedly
- **CryLasLaserError** – if laser\_off() or close\_shutter() fail

**property target\_pulse\_energy**

**update\_laser\_status()** → *None*

Update the laser status to *LaserStatus.NOT\_READY* or *LaserStatus.INACTIVE* or *LaserStatus.ACTIVE*.

Note: laser never explicitly says that it is not ready (*LaserStatus.NOT\_READY*) in response to ‘STATUS’ command. It only says that it is ready (heated-up and implicitly inactive/off) or active (on). If it’s not either of these then the answer is *Answers.HEAD*. Moreover, the only time the laser explicitly says that its status is inactive (*Answers.INACTIVE*) is after issuing a ‘LASER OFF’ command.

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened

**update\_repetition\_rate()** → *None*

Query the laser repetition rate.

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

**update\_shutter\_status()** → *None*

Update the shutter status (OPENED or CLOSED)

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

**update\_target\_pulse\_energy()** → *None*

Query the laser pulse energy.

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

**wait\_until\_ready()** → *None*

Block execution until the laser is ready

**Raises**

- **CryLasLaserError** – if the polling thread stops before the laser is ready

```
class CryLasLaserConfig(calibration_factor: Union[int, float] = 4.35, polling_period: Union[int, float] = 12,
                        polling_timeout: Union[int, float] = 300, auto_laser_on: bool = True,
                        init_shutter_status: Union[int, CryLasLaserShutterStatus] =
                        CryLasLaserShutterStatus.CLOSED)
```

Bases: `object`

Device configuration dataclass for the CryLas laser controller.

#### **ShutterStatus**

alias of `CryLasLaserShutterStatus`

**auto\_laser\_on:** `bool = True`

**calibration\_factor:** `Union[int, float] = 4.35`

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### **Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**init\_shutter\_status:** `Union[int, CryLasLaserShutterStatus] = 0`

**is\_configdataclass** = `True`

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

#### **Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### **Returns**

a list of strings containing all optional keys.

**polling\_period:** `Union[int, float] = 12`

**polling\_timeout:** `Union[int, float] = 300`

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

#### **Returns**

a list of strings containing all required keys.

**exception CryLasLaserError**

Bases: `DeviceError`

General error with the CryLas Laser.

**exception CryLasLaserNotReadyError**Bases: *CryLasLaserError*

Error when trying to turn on the CryLas Laser before it is ready.

**class CryLasLaserPoller(spoll\_handler: Callable, check\_handler: Callable, check\_laser\_status\_handler: Callable, polling\_delay\_sec: Union[int, float] = 0, polling\_interval\_sec: Union[int, float] = 1, polling\_timeout\_sec: Optional[Union[int, float]] = None)**Bases: *Poller*

Poller class for polling the laser status until the laser is ready.

**Raises**

- *CryLasLaserError* – if the timeout is reached before the laser is ready
- *SerialCommunicationIOError* – when communication port is closed.

**class CryLasLaserSerialCommunication(configuration)**Bases: *SerialCommunication*

Specific communication protocol implementation for the CryLas laser controller. Already predefines device-specific protocol parameters in config.

**READ\_TEXT\_SKIP\_PREFIXES = ('>', 'MODE:')**

Prefixes of lines that are skipped when read from the serial port.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**query(cmd: str, prefix: str, post\_cmd: Optional[str] = None) → str**

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns the line in question.

**Parameters**

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer
- **post\_cmd** – optional additional command to send after the query

**Returns**

line in question as a string

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**query\_all(cmd: str, prefix: str)**

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns a list of successive lines starting with prefix.

**Parameters**

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer

**Returns**

line in question as a string

**Raises**

`SerialCommunicationIOError` – when communication port is not opened

**read()** → str

Read first line of text from the serial port that does not start with any of `self.READ_TEXT_SKIP_PREFIXES`.

**Returns**

String read from the serial port; '' if there was nothing to read.

**Raises**

`SerialCommunicationIOError` – when communication port is not opened

```
class CryLasLaserSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Union[str, NoneType] = None, baudrate: int = 19200,
                                             parity: Union[str,
                                             hvl_ccb.comm.serial.SerialCommunicationParity] =
                                             <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                             Union[int,
                                             hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                             <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                             Union[int,
                                             hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                             <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                             Union[int, float] = 10)
```

Bases: `SerialCommunicationConfig`

**baudrate: int = 19200**

Baudrate for CryLas laser is 19200 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

```

parity: Union[str, SerialCommunicationParity] = 'N'
    CryLas laser does not use parity

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns
    a list of strings containing all required keys.

stopbits: Union[int, SerialCommunicationStopbits] = 1
    CryLas laser uses one stop bit

terminator: bytes = b'\n'
    The terminator is LF

timeout: Union[int, float] = 10
    use 10 seconds timeout as default (a long timeout is needed!)


class CryLasLaserShutterStatus(value=<no_arg>, names=None, module=None, qualname=None,
    type=None, start=1, boundary=None)

Bases: Enum
Status of the CryLas laser shutter

CLOSED = 0
OPENED = 1

```

## [hvl\\_ccb.dev.ea\\_psi9000](#)

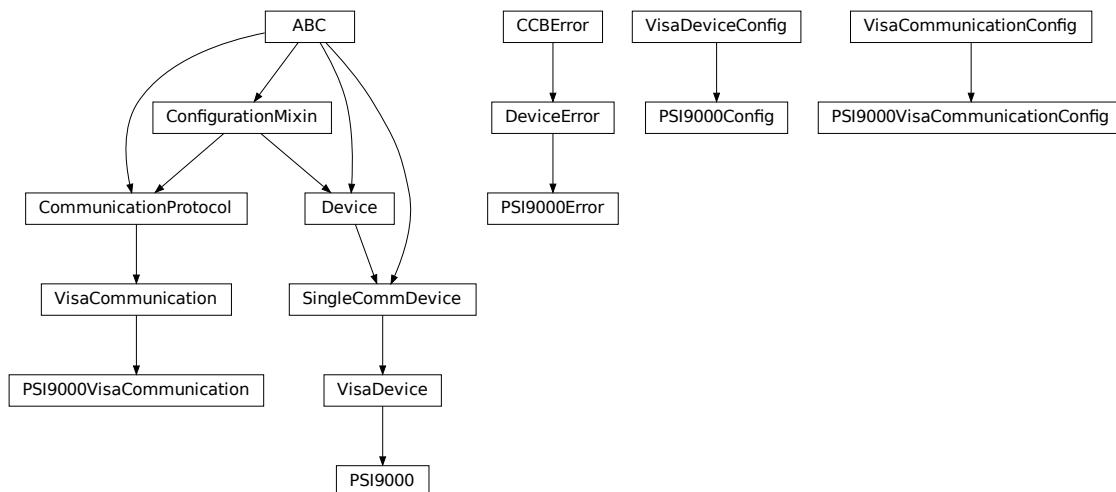

Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to now, all the testing was done with NI-Visa)

```
class PSI9000(com: Union[PSI9000VisaCommunication, PSI9000VisaCommunicationConfig, dict], dev_config:  
    Optional[Union[PSI9000Config, dict]] = None)
```

Bases: *VisaDevice*

Elektro Automatik PSI 9000 power supply.

**MS\_NOMINAL\_CURRENT = 2040**

**MS\_NOMINAL\_VOLTAGE = 80**

**SHUTDOWN\_CURRENT\_LIMIT = 0.1**

**SHUTDOWN\_VOLTAGE\_LIMIT = 0.1**

**check\_master\_slave\_config()** → *None*

Checks if the master / slave configuration and initializes if successful

**Returns**

*PSI9000Error* – if master-slave configuration failed

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls()**

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns**

the VisaCommunication class

**get\_output()** → bool

Reads the current state of the DC output of the source. Returns True, if it is enabled, false otherwise.

**Returns**

the state of the DC output

**get\_system\_lock()** → bool

Get the current lock state of the system. The lock state is true, if the remote control is active and false, if not.

**Returns**

the current lock state of the device

**get\_ui\_lower\_limits()** → Tuple[float, float]

Get the lower voltage and current limits. A lower power limit does not exist.

**Returns**

Umin in V, Imin in A

**get\_uip\_upper\_limits()** → Tuple[float, float, float]

Get the upper voltage, current and power limits.

**Returns**

Umax in V, Imax in A, Pmax in W

**get\_voltage\_current\_setpoint()** → Tuple[float, float]

Get the voltage and current setpoint of the current source.

**Returns**

Uset in V, Iset in A

**measure\_voltage\_current()** → Tuple[float, float]

Measure the DC output voltage and current

**Returns**

Umeas in V, Imeas in A

**set\_lower\_limits(voltage\_limit: Optional[float] = None, current\_limit: Optional[float] = None)** → *None*

Set the lower limits for voltage and current. After writing the values a check is performed if the values are set correctly.

**Parameters**

- **voltage\_limit** – is the lower voltage limit in V
- **current\_limit** – is the lower current limit in A

**Raises**

**PSI9000Error** – if the limits are out of range

**set\_output(target\_onstate: bool)** → *None*

Enables / disables the DC output.

**Parameters**

**target\_onstate** – enable or disable the output power

**Raises**

**PSI9000Error** – if operation was not successful

**set\_system\_lock(lock: bool)** → *None*

Lock / unlock the device, after locking the control is limited to this class unlocking only possible when voltage and current are below the defined limits

**Parameters**

**lock** – True: locking, False: unlocking

**set\_upper\_limits(voltage\_limit: Optional[float] = None, current\_limit: Optional[float] = None, power\_limit: Optional[float] = None)** → *None*

Set the upper limits for voltage, current and power. After writing the values a check is performed if the values are set. If a parameter is left blank, the maximum configurable limit is set.

**Parameters**

- **voltage\_limit** – is the voltage limit in V
- **current\_limit** – is the current limit in A
- **power\_limit** – is the power limit in W

**Raises**

**PSI9000Error** – if limits are out of range

**set\_voltage\_current(volt: float, current: float)** → *None*

Set voltage and current setpoints.

After setting voltage and current, a check is performed if writing was successful.

**Parameters**

- **volt** – is the setpoint voltage: 0.81.6 V (1.02 \* 0-80 V) (absolute max, can be smaller if limits are set)
- **current** – is the setpoint current: 0..2080.8 A (1.02 \* 0 - 2040 A) (absolute max, can be smaller if limits are set)

**Raises**

**PSI9000Error** – if the desired setpoint is out of limits

**start()** → *None*

Start this device.

**stop()** → *None*

Stop this device. Turns off output and lock, if enabled.

```
class PSI9000Config(spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2,
                     power_limit: Union[int, float] = 43500, voltage_lower_limit: Union[int, float] = 0.0,
                     voltage_upper_limit: Union[int, float] = 10.0, current_lower_limit: Union[int, float] = 0.0,
                     current_upper_limit: Union[int, float] = 2040.0, wait_sec_system_lock: Union[int, float]
                     = 0.5, wait_sec_settings_effect: Union[int, float] = 1, wait_sec_initialisation: Union[int,
                     float] = 2)
```

Bases: *VisaDeviceConfig*

Elektro Automatik PSI 9000 power supply device class. The device is communicating over a VISA TCP socket.

Using this power supply, DC voltage and current can be supplied to a load with up to 2040 A and 80 V (using all four available units in parallel). The maximum power is limited by the grid, being at 43.5 kW available through the CEE63 power socket.

**clean\_values()** → *None*

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**current\_lower\_limit: Union[int, float] = 0.0**

Lower current limit in A, depending on the experimental setup.

**current\_upper\_limit: Union[int, float] = 2040.0**

Upper current limit in A, depending on the experimental setup.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**power\_limit: Union[int, float] = 43500**

Power limit in W depending on the experimental setup. With 3x63A, this is 43.5kW. Do not change this value, if you do not know what you are doing. There is no lower power limit.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**voltage\_lower\_limit: Union[int, float] = 0.0**

Lower voltage limit in V, depending on the experimental setup.

**voltage\_upper\_limit: Union[int, float] = 10.0**

Upper voltage limit in V, depending on the experimental setup.

**wait\_sec\_initialisation: Union[int, float] = 2****wait\_sec\_settings\_effect: Union[int, float] = 1****wait\_sec\_system\_lock: Union[int, float] = 0.5****exception PSI9000Error**

Bases: *DeviceError*

Base error class regarding problems with the PSI 9000 supply.

**class PSI9000VisaCommunication(*configuration*)**

Bases: *VisaCommunication*

Communication protocol used with the PSI 9000 power supply.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class PSI9000VisaCommunicationConfig(*host: Union[str, IPv4Address, IPv6Address], interface\_type: Union[str, InterfaceType] = InterfaceType.TCPIP\_SOCKET, board: int = 0, port: int = 5025, timeout: int = 5000, chunk\_size: int = 204800, open\_timeout: int = 1000, write\_termination: str = '\n', read\_termination: str = '\n', visa\_backend: str = ''*)**

Bases: *VisaCommunicationConfig*

Visa communication protocol config dataclass with specification for the PSI 9000 power supply.

**force\_value(*fieldname, value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**interface\_type:** Union[str, *InterfaceType*] = 1

Interface type of the VISA connection, being one of *InterfaceType*.

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.


**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

## hvl\_ccb.dev.fug

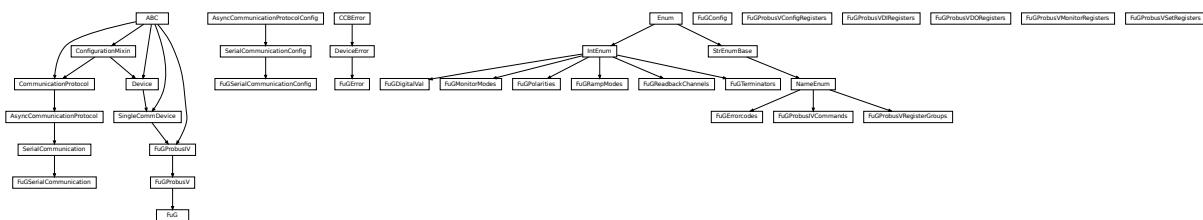

Device classes for “Probus V - ADDAT30” Interfaces which are used to control power supplies from FuG Elektronik GmbH

This interface is used for many FuG power units. Manufacturer homepage: <https://www.fug-elektronik.de>

The Professional Series of Power Supplies from FuG is a series of low, medium and high voltage direct current power supplies as well as capacitor chargers. The class FuG is tested with a HCK 800-20 000 in Standard Mode. The addressable mode is not implemented. Check the code carefully before using it with other devices. Manufacturer homepage: <https://www.fug-elektronik.de/netzgeraete/professional-series/>

The documentation of the interface from the manufacturer can be found here: [https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus\\_V.zip](https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus_V.zip)

The provided classes support the basic and some advanced commands. The commands for calibrating the power supplies are not implemented, as they are only for very special porpoises and should not be used by “normal” customers.

**class FuG(com, dev\_config=None)**

Bases: *FuGProbusV*

FuG power supply device class.

The power supply is controlled over a FuG ADDA Interface with the PROBUS V protocol

**property config\_status:** *FuGProbusVConfigRegisters*

Returns the registers for the registers with the configuration and status values

**Returns**

FuGProbusVConfigRegisters

**property current:** *FuGProbusVSetRegisters*

Returns the registers for the current output

**Returns**

**property current\_monitor:** *FuGProbusVMonitorRegisters*

Returns the registers for the current monitor.

A typically usage will be “self.current\_monitor.value” to measure the output current

**Returns**

**property di:** *FuGProbusVDIRegisters*

Returns the registers for the digital inputs

**Returns**

FuGProbusVDIRegisters

**identify\_device()** → *None*

Identify the device nominal voltage and current based on its model number.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**property max\_current:** *Union[int, float]*

Returns the maximal current which could provided within the test setup

**Returns**

**property max\_current\_hardware:** *Union[int, float]*

Returns the maximal current which could provided with the power supply

**Returns**

**property max\_voltage:** *Union[int, float]*

Returns the maximal voltage which could provided within the test setup

**Returns**

**property max\_voltage\_hardware:** *Union[int, float]*

Returns the maximal voltage which could provided with the power supply

**Returns**

**property on:** *FuGProbusVDORegisters*

Returns the registers for the output switch to turn the output on or off

**Returns**

FuGProbusVDORegisters

**property outX0:** *FuGProbusVDORegisters*

Returns the registers for the digital output X0

**Returns**

FuGProbusVDORegisters

**property outX1: *FuGProbusVDORegisters***

Returns the registers for the digital output X1

**Returns**

FuGProbusVDORegisters

**property outX2: *FuGProbusVDORegisters***

Returns the registers for the digital output X2

**Returns**

FuGProbusVDORegisters

**property outXCMD: *FuGProbusVDORegisters***

Returns the registers for the digital outputX-CMD

**Returns**

FuGProbusVDORegisters

**start(*max\_voltage=0, max\_current=0*) → *None***

Opens the communication protocol and configures the device.

**Parameters**

- **max\_voltage** – Configure here the maximal permissible voltage which is allowed in the given experimental setup
- **max\_current** – Configure here the maximal permissible current which is allowed in the given experimental setup

**property voltage: *FuGProbusVSetRegisters***

Returns the registers for the voltage output

**Returns**

**property voltage\_monitor: *FuGProbusVMonitorRegisters***

Returns the registers for the voltage monitor.

A typically usage will be “self.voltage\_monitor.value” to measure the output voltage

**Returns**

**class FuGConfig(*wait\_sec\_stop\_commands: Union[int, float] = 0.5*)**

Bases: object

Device configuration dataclass for FuG power supplies.

**clean\_values()**

**force\_value(*fieldname, value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

```
classmethod optional_defaults() → Dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

```
wait_sec_stop_commands: Union[int, float] = 0.5
```

Time to wait after subsequent commands during stop (in seconds)

```
class FuGDigitalVal(value)
```

Bases: IntEnum

An enumeration.

```
NO = 0
```

```
OFF = 0
```

```
ON = 1
```

```
YES = 1
```

```
exception FuGError(*args, **kwargs)
```

Bases: *DeviceError*

Error with the FuG voltage source.

```
errorcode: str
```

Errorcode from the Probus, see documentation of Probus V chapter 5. Errors with three-digit errorcodes are thrown by this python module.

```
class FuGErrorcodes(value=<no_arg>, name=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: *NameEnum*

The power supply can return an errorcode. These errorcodes are handled by this class. The original errorcodes from the source are with one or two digits, see documentation of Probus V chapter 5. All three-digit errorcodes are from this python module.

```
E0 = ('no error', 'standard response on each command')
```

```
E1 = ('no data available', 'Customer tried to read from GPIB but there were no data prepared. (IBIG50 sent command ~T2 to ADDA)')
```

```
E10 = ('unknown SCPI command', 'This SCPI command is not implemented')
```

```
E100 = ('Command is not implemented', 'You tried to execute a command, which is not implemented or does not exist')

E106 = ('The rampstate is a read-only register', 'You tried to write data to the register, which can only give you the status of the ramping.')

E11 = ('not allowed Trigger-on-Talk', 'Not allowed attempt to Trigger-on-Talk (~T1) while ADDA was in addressable mode.')

E115 = ('The given index to select a digital value is out of range', 'Only integer values between 0 and 1 are allowed.')

E12 = ('invalid argument in ~Tn command', 'Only ~T1 and ~T2 is implemented.')

E125 = ('The given index to select a ramp mode is out of range', 'Only integer values between 0 and 4 are allowed.')

E13 = ('invalid N-value', 'Register > K8 contained an invalid value. Error code is output on an attempt to query data with ? or ~T1')

E135 = ('The given index to select the readback channel is out of range', 'Only integer values between 0 and 6 are allowed.')

E14 = ('register is write only', 'Some registers can only be written to (i.e.> H0)')

E145 = ('The given value for the AD-conversion is unknown', 'Valid values for the ad-conversion are integer values from "0" to "7".')

E15 = ('string too long', 'i.e. serial number string too long during calibration')

E155 = ('The given value to select a polarity is out range.', 'The value should be 0 or 1.')

E16 = ('wrong checksum', 'checksum over command string was not correct, refer also to 4.4 of the Probus V documentation')

E165 = ('The given index to select the terminator string is out of range', '')

E2 = ('unknown register type', "No valid register type after '>'")

E206 = ('This status register is read-only', 'You tried to write data to this register, which can only give you the actual status of the corresponding digital output.')

E306 = ('The monitor register is read-only', 'You tried to write data to a monitor, which can only give you measured data.')

E4 = ('invalid argument', 'The argument of the command was rejected .i.e. malformed number')

E5 = ('argument out of range', 'i.e. setvalue higher than type value')

E504 = ('Empty string as response', 'The connection is broken.')

E505 = ('The returned register is not the requested.', 'Maybe the connection is overburden.')
```

```
E6 = ('register is read only', 'Some registers can only be read but not written to.  
(i.e. monitor registers)')  
  
E666 = ('You cannot overwrite the most recent error in the interface of the power  
supply. But, well: You created an error anyway...', '')  
  
E7 = ('Receive Overflow', 'Command string was longer than 50 characters.')  
  
E8 = ('EEPROM is write protected', 'Write attempt to calibration data while the  
write protection switch was set to write protected.')  
  
E9 = ('address error', 'A non addressed command was sent to ADDA while it was in  
addressable mode (and vice versa).')  
  
raise_()  
  
class FuGMonitorModes(value)  
    Bases: IntEnum  
    An enumeration.  
  
    T1MS = 1  
        15 bit + sign, 1 ms integration time  
    T200MS = 6  
        typ. 19 bit + sign, 200 ms integration time  
    T20MS = 3  
        17 bit + sign, 20 ms integration time  
    T256US = 0  
        14 bit + sign, 256 us integration time  
    T40MS = 4  
        17 bit + sign, 40 ms integration time  
    T4MS = 2  
        15 bit + sign, 4 ms integration time  
    T800MS = 7  
        typ. 20 bit + sign, 800 ms integration time  
    T80MS = 5  
        typ. 18 bit + sign, 80 ms integration time  
  
class FuGPolarities(value)  
    Bases: IntEnum  
    An enumeration.  
    NEGATIVE = 1  
    POSITIVE = 0  
  
class FuGProbusIV(com, dev_config=None)  
    Bases: SingleCommDevice, ABC  
    FuG Probus IV device class  
    Sends basic SCPI commands and reads the answer. Only the special commands and PROBUS IV instruction set  
    is implemented.
```

**command**(*command*: FuGProbusIVCommands, *value*=None) → str

**Returns**

- **command** – one of the commands given within FuGProbusIVCommands
- **value** – an optional value, depending on the command

**Returns**

a String if a query was performed

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**output\_off()** → *None*

Switch DC voltage output off.

**reset()** → *None*

Reset of the interface: All setvalues are set to zero

**abstract start()**

Open the associated communication protocol.

**stop()** → *None*

Close the associated communication protocol.

**class FuGProbusIVCommands**(*value*=<no\_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *NameEnum*

An enumeration.

**ADMODE** = ('S', (<enum 'FuGMonitorModes'>, <class 'int'>))

**CURRENT** = ('I', (<class 'int'>, <class 'float'>))

**EXECUTE** = ('X', None)

**EXECUTEONX** = ('G', (<enum 'FuGDigitalVal'>, <class 'int'>))

Wait for "X" to execute pending commands

**ID** = ('\*IDN?', None)

**OUTPUT** = ('F', (<enum 'FuGDigitalVal'>, <class 'int'>))

**POLARITY** = ('P', (<enum 'FuGPolarities'>, <class 'int'>))

**QUERY** = ('?', None)

**READBACKCHANNEL** = ('N', (<enum 'FuGReadbackChannels'>, <class 'int'>))

**RESET** = ('=', None)

```

TERMINATOR = ('Y', (<enum 'FuGTerminators'>, <class 'int'>))
VOLTAGE = ('U', (<class 'int'>, <class 'float'>))
XOUTPUTS = ('R', <class 'int'>)
TODO: the possible values are limited to 0..13

class FuGProbusV(com, dev_config=None)
Bases: FuGProbusIV
FuG Probus V class which uses register based commands to control the power supplies
get_register(register: str) → str
    get the value from a register
Parameters
    register – the register from which the value is requested
Returns
    the value of the register as a String
set_register(register: str, value: Union[int, float, str]) → None
    generic method to set value to register
Parameters
    • register – the name of the register to set the value
    • value – which should be written to the register

class FuGProbusVConfigRegisters(fug, super_register: FuGProbusVRegisterGroups)
Bases: object
Configuration and Status values, acc. 4.2.5
property execute_on_x: FuGDigitalVal
    status of Execute-on-X
Returns
    FuGDigitalVal of the status
property most_recent_error: FuGErrorcodes
    Reads the Error-Code of the most recent command
Return FuGError
Raises
    FuGError – if code is not “E0”
property readback_data: FuGReadbackChannels
    Preselection of readout data for Trigger-on-Talk
Returns
    index for the readback channel
property srq_mask: int
    SRQ-Mask, Service-Request Enable status bits for SRQ 0: no SRQ Bit 2: SRQ on change of status to CC
    Bit 1: SRQ on change to CV
Returns
    representative integer value

```

```
property srq_status: str
    SRQ-Statusbyte output as a decimal number: Bit 2: PS is in CC mode Bit 1: PS is in CV mode

    Returns
        representative string

property status: str
    Statusbyte as a string of 0/1. Combined status (compatibel to Probus IV), MSB first: Bit 7: I-REG Bit 6: V-REG Bit 5: ON-Status Bit 4: 3-Reg Bit 3: X-Stat (polarity) Bit 2: Cal-Mode Bit 1: unused Bit 0: SEL-D

    Returns
        string of 0/1

property terminator: FuGTerminators
    Terminator character for answer strings from ADDA

    Returns
        FuGTerminators

class FuGProbusVDIRegisters(fug, super_register: FuGProbusVRegisterGroups)
    Bases: object

    Digital Inputs acc. 4.2.4

    property analog_control: FuGDigitalVal

        Returns
            shows 1 if power supply is controlled by the analog interface

    property calibration_mode: FuGDigitalVal

        Returns
            shows 1 if power supply is in calibration mode

    property cc_mode: FuGDigitalVal

        Returns
            shows 1 if power supply is in CC mode

    property cv_mode: FuGDigitalVal

        Returns
            shows 1 if power supply is in CV mode

    property digital_control: FuGDigitalVal

        Returns
            shows 1 if power supply is digitally controlled

    property on: FuGDigitalVal

        Returns
            shows 1 if power supply ON

    property reg_3: FuGDigitalVal

        For special applications.

        Returns
            input from bit 3-REG
```

```
property x_stat: FuGPolarities
```

**Returns**

polarity of HVPS with polarity reversal

```
class FuGProbusVDORegisters(fug, super_register: FuGProbusVRegisterGroups)
```

Bases: object

Digital outputs acc. 4.2.2

```
property out: Union[int, FuGDigitalVal]
```

Status of the output according to the last setting. This can differ from the actual state if output should only pulse.

**Returns**

FuGDigitalVal

```
property status: FuGDigitalVal
```

Returns the actual value of output. This can differ from the set value if pulse function is used.

**Returns**

FuGDigitalVal

```
class FuGProbusVMonitorRegisters(fug, super_register: FuGProbusVRegisterGroups)
```

Bases: object

Analog monitors acc. 4.2.3

```
property adc_mode: FuGMonitorModes
```

The programmed resolution and integration time of the AD converter

**Returns**

FuGMonitorModes

```
property value: float
```

Value from the monitor.

**Returns**

a float value in V or A

```
property value_raw: float
```

uncalibrated raw value from AD converter

**Returns**

float value from ADC

```
class FuGProbusVRegisterGroups(value=<no_arg>, names=None, module=None, qualname=None,  
type=None, start=1, boundary=None)
```

Bases: *NameEnum*

An enumeration.

**CONFIG = 'K'**

**INPUT = 'D'**

**MONITOR\_I = 'M1'**

**MONITOR\_V = 'M0'**

**OUTPUTONCMD = 'BON'**

```
OUTPUTX0 = 'B0'
OUTPUTX1 = 'B1'
OUTPUTX2 = 'B2'
OUTPUTXCMD = 'BX'
SETCURRENT = 'S1'
SETVOLTAGE = 'S0'

class FuGProbusVSetRegisters(fug, super_register: FuGProbusVRegisterGroups)
Bases: object

Setvalue control acc. 4.2.1 for the voltage and the current output

property actualsetvalue: float
    The actual valid set value, which depends on the ramp function.

    Returns
        actual valid set value

property high_resolution: FuGDigitalVal
    Status of the high resolution mode of the output.

    Return 0
        normal operation

    Return 1
        High Res. Mode

property rampmode: FuGRampModes
    The set ramp mode to control the setvalue.

    Returns
        the mode of the ramp as instance of FuGRampModes

property ramprate: float
    The set ramp rate in V/s.

    Returns
        ramp rate in V/s

property rampstate: FuGDigitalVal
    Status of ramp function.

    Return 0
        if final setvalue is reached

    Return 1
        if still ramping up

property setvalue: float
    For the voltage or current output this setvalue was programmed.

    Returns
        the programmed setvalue
```

```
class FuGRampModes(value)
```

Bases: IntEnum

An enumeration.

**FOLLOWRAMP = 1**

Follow the ramp up- and downwards

**IMMEDIATELY = 0**

Standard mode: no ramp

**ONLYUPWARDSOFFTOZERO = 4**

Follow the ramp up- and downwards, if output is OFF set value is zero

**RAMPUPWARDS = 2**

Follow the ramp only upwards, downwards immediately

**SPECIALRAMPUPWARDS = 3**

Follow a special ramp function only upwards

```
class FuGReadbackChannels(value)
```

Bases: IntEnum

An enumeration.

**CURRENT = 1**

**FIRMWARE = 5**

**RATEDCURRENT = 4**

**RATEDVOLTAGE = 3**

**SN = 6**

**STATUSBYTE = 2**

**VOLTAGE = 0**

```
class FuGSerialCommunication(configuration)
```

Bases: *SerialCommunication*

Specific communication protocol implementation for FuG power supplies. Already defines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**query(*command: str*) → str**

Send a command to the interface and handle the status message. Raises an error, if the answer starts with “E”.

**Parameters**

**command** – Command to send

**Raises**

**FuError** – if the connection is broken or the error from the power source itself

**Returns**

Answer from the interface or empty string

```
class FuGSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                    encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                    Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int =
                                    10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
                                    <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
                                    hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                    <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
                                    hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int,
                                    float] = 3)
```

Bases: *SerialCommunicationConfig*

**baudrate: int = 9600**

Baudrate for FuG power supplies is 9600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**default\_n\_attempts\_read\_text\_nonempty: int = 10**

default number of attempts to read a non-empty text

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

FuG does not use parity

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.


```

stopbits: Union[int, SerialCommunicationStopbits] = 1
    FuG uses one stop bit
terminator: bytes = b'\n'
    The terminator is LF
timeout: Union[int, float] = 3
    use 3 seconds timeout as default
wait_sec_read_text_nonempty: Union[int, float] = 0.5
    default time to wait between attempts of reading a non-empty text

class FuGTerminators(value)
    Bases: IntEnum
    An enumeration.
    CR = 3
    CRLF = 0
    LF = 2
    LFCR = 1

```

## hvl\_ccb.dev.heinzinger

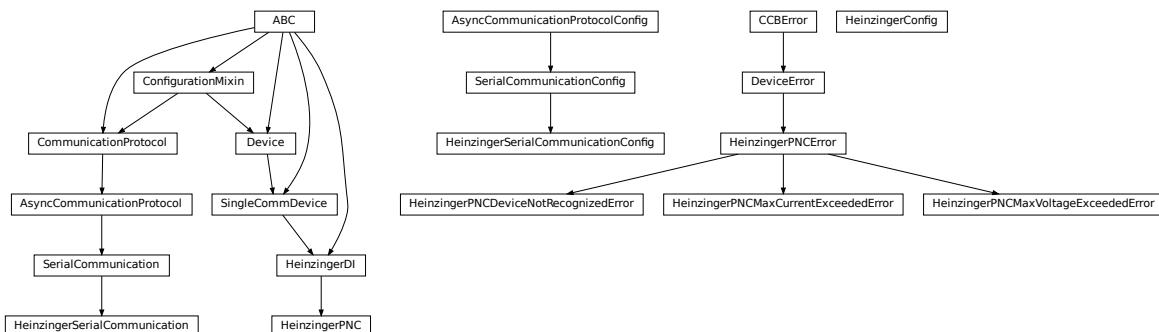

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Manufacturer homepage: <https://www.heinzinger.com/products/accessories-and-more/digital-interfaces/>

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class **HeinzingerPNC** is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCCap. Manufacturer homepage: <https://www.heinzinger.com/products/high-voltage/universal-high-voltage-power-supplies/>

```

class HeinzingerConfig(default_number_of_recordings: Union[int, RecordingsEnum] = 1,
number_of_decimals: int = 6, wait_sec_stop_commands: Union[int, float] = 0.5)

Bases: object
Device configuration dataclass for Heinzinger power supplies.

```

```
class RecordingsEnum(value)
    Bases: IntEnum
    An enumeration.

    EIGHT = 8
    FOUR = 4
    ONE = 1
    SIXTEEN = 16
    TWO = 2

    clean_values()

    default_number_of_recordings: Union[int, RecordingsEnum] = 1

    force_value(fieldname, value)
        Forces a value to a dataclass field despite the class being frozen.

        NOTE: you can define post_force_value method with same signature as this method to do extra processing
        after value has been forced on fieldname.

    Parameters
        • fieldname – name of the field
        • value – value to assign

    is_configdataclass = True

    classmethod keys() → Sequence[str]
        Returns a list of all configdataclass fields key-names.

    Returns
        a list of strings containing all keys.

    number_of_decimals: int = 6

    classmethod optional_defaults() → Dict[str, object]
        Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
        on instantiation.

    Returns
        a list of strings containing all optional keys.

    classmethod required_keys() → Sequence[str]
        Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
        instantiation.

    Returns
        a list of strings containing all required keys.

    wait_sec_stop_commands: Union[int, float] = 0.5
        Time to wait after subsequent commands during stop (in seconds)
```

```
class HeinzingerDI(com, dev_config=None)
Bases: SingleCommDevice, ABC
Heinzinger Digital Interface I/II device class
Sends basic SCPI commands and reads the answer. Only the standard instruction set from the manual is implemented.

class OutputStatus(value)
    Bases: IntEnum
    Status of the voltage output
    OFF = 0
    ON = 1
    UNKNOWN = -1

static config_cls()
    Return the default configdataclass class.

    Returns
        a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns
        the type of the standard communication protocol for this device

get_current() → float
    Queries the set current of the Heinzinger PNC (not the measured current!).

    Raises
        SerialCommunicationIOError – when communication port is not opened

get_interface_version() → str
    Queries the version number of the digital interface.

    Raises
        SerialCommunicationIOError – when communication port is not opened

get_number_of_recordings() → int
    Queries the number of recordings the device is using for average value calculation.

    Returns
        int number of recordings

    Raises
        SerialCommunicationIOError – when communication port is not opened

get_serial_number() → str
    Ask the device for its serial number and returns the answer as a string.

    Returns
        string containing the device serial number

    Raises
        SerialCommunicationIOError – when communication port is not opened
```

**get\_voltage()** → float

Queries the set voltage of the Heinzinger PNC (not the measured voltage!).

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**measure\_current()** → float

Ask the Device to measure its output current and return the measurement result.

**Returns**

measured current as float

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**measure\_voltage()** → float

Ask the Device to measure its output voltage and return the measurement result.

**Returns**

measured voltage as float

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**output\_off()** → *None*

Switch DC voltage output off and updates the output status.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**output\_on()** → *None*

Switch DC voltage output on and updates the output status.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**property output\_status: OutputStatus**

**reset\_interface()** → *None*

Reset of the digital interface; only Digital Interface I: Power supply is switched to the Local-Mode (Manual operation)

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**set\_current(value: Union[int, float])** → *None*

Sets the output current of the Heinzinger PNC to the given value.

**Parameters**

**value** – current expressed in *self.unit\_current*

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**set\_number\_of\_recordings(value: Union[int, RecordingsEnum])** → *None*

Sets the number of recordings the device is using for average value calculation. The possible values are 1, 2, 4, 8 and 16.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**set\_voltage**(*value*: Union[int, float]) → *None*

Sets the output voltage of the Heinzinger PNC to the given value.

**Parameters**

**value** – voltage expressed in *self.unit\_voltage*

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**abstract start()**

Opens the communication protocol.

**Raises**

*SerialCommunicationIOError* – when communication port cannot be opened.

**stop()** → *None*

Stop the device. Closes also the communication protocol.

**class HeinzingerPNC**(*com*, *dev\_config*=None)

Bases: *HeinzingerDI*

Heinzinger PNC power supply device class.

The power supply is controlled over a Heinzinger Digital Interface I/II

**class UnitCurrent**(*value*=<no\_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *AutoNumberNameEnum*

An enumeration.

**A** = 3

**UNKNOWN** = 1

**mA** = 2

**class UnitVoltage**(*value*=<no\_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *AutoNumberNameEnum*

An enumeration.

**UNKNOWN** = 1

**V** = 2

**kV** = 3

**identify\_device()** → *None*

Identify the device nominal voltage and current based on its serial number.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**property max\_current:** Union[int, float]

**property max\_current\_hardware:** Union[int, float]

**property max\_voltage:** Union[int, float]

**property max\_voltage\_hardware:** Union[int, float]

**set\_current**(value: Union[int, float]) → *None*

Sets the output current of the Heinzinger PNC to the given value.

**Parameters**

**value** – current expressed in *self.unit\_current*

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**set\_voltage**(value: Union[int, float]) → *None*

Sets the output voltage of the Heinzinger PNC to the given value.

**Parameters**

**value** – voltage expressed in *self.unit\_voltage*

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**start()** → *None*

Opens the communication protocol and configures the device.

**property unit\_current:** UnitCurrent

**property unit\_voltage:** UnitVoltage

**exception HeinzingerPNCDeviceNotRecognizedError**

Bases: *HeinzingerPNCError*

Error indicating that the serial number of the device is not recognized.

**exception HeinzingerPNCError**

Bases: *DeviceError*

General error with the Heinzinger PNC voltage source.

**exception HeinzingerPNCMaxCurrentExceededError**

Bases: *HeinzingerPNCError*

Error indicating that program attempted to set the current to a value exceeding ‘max\_current’.

**exception HeinzingerPNCMaxVoltageExceededError**

Bases: *HeinzingerPNCError*

Error indicating that program attempted to set the voltage to a value exceeding ‘max\_voltage’.

**class HeinzingerSerialCommunication**(*configuration*)

Bases: *SerialCommunication*

Specific communication protocol implementation for Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
class HeinzingerSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
    encoding_error_handling: str = 'strict',
    wait_sec_read_text_nonempty: Union[int, float] = 0.5,
    default_n_attempts_read_text_nonempty: int = 40, port:
    Union[str, NoneType] = None, baudrate: int = 9600, parity:
    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
    <SerialCommunicationParity.NONE: 'N'>, stopbits:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationStopbits] =
    <SerialCommunicationStopbits.ONE: 1>, bytesize:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationBytesize] =
    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
    Union[int, float] = 3)
```

Bases: *SerialCommunicationConfig*

**baudrate: int = 9600**

Baudrate for Heinzinger power supplies is 9600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**default\_n\_attempts\_read\_text\_nonempty: int = 40**

increased to 40 default number of attempts to read a non-empty text

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

#### Returns

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### Returns

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

Heinzinger does not use parity

**classmethod required\_keys() → Sequence[str]**


Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

#### Returns

a list of strings containing all required keys.

```
stopbits: Union[int, SerialCommunicationStopbits] = 1
    Heinzinger uses one stop bit
terminator: bytes = b'\n'
    The terminator is LF
timeout: Union[int, float] = 3
    use 3 seconds timeout as default
wait_sec_read_text_nonempty: Union[int, float] = 0.5
    default time to wait between attempts of reading a non-empty text
```

## hvl\_ccb.dev.labjack

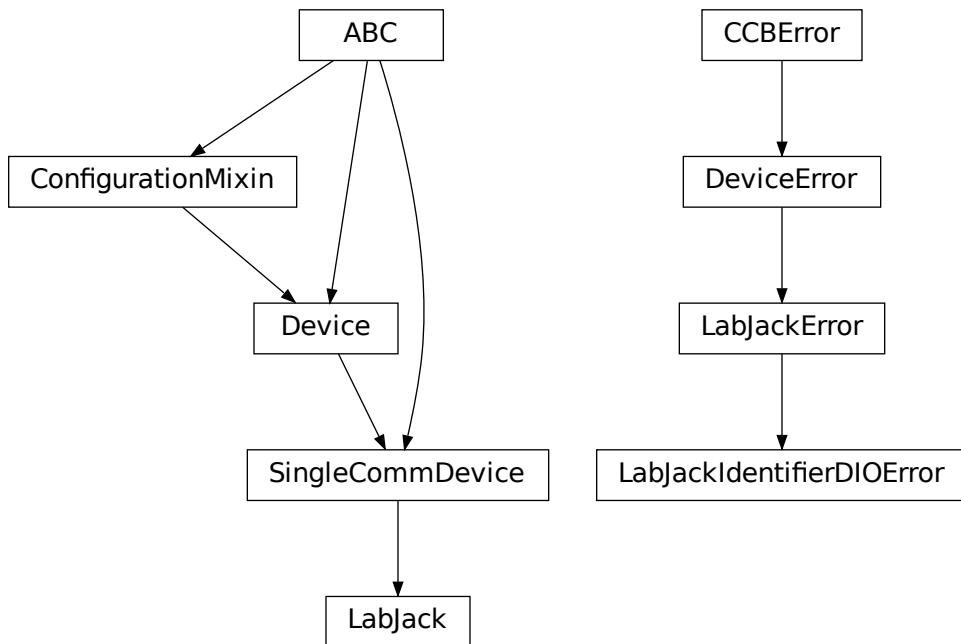

A  
LabJack T-series devices wrapper around the LabJack's LJM Library; see <https://labjack.com/ljm> . The wrapper was originally developed and tested for a LabJack T7-PRO device.

## Extra installation

To use this LabJack T-series devices wrapper:

1. install the hvl\_ccb package with a labjack extra feature:

```
$ pip install "hvl_ccb[labjack]"
```

this will install the Python bindings for the library.

2. install the library - follow instruction in <https://labjack.com/support/software/installers/ljm> .

**class LabJack**(*com*, *dev\_config=None*)

Bases: *SingleCommDevice*

LabJack Device.

This class is tested with a LabJack T7-Pro and should also work with T4 and T7 devices communicating through the LJM Library. Other or older hardware versions and variants of LabJack devices are not supported.

**class AInRange**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *StrEnumBase*

An enumeration.

**ONE = 1.0**

**ONE\_HUNDREDTH = 0.01**

**ONE\_TENTH = 0.1**

**TEN = 10.0**

**property value: float**

**class BitLimit**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *IntEnum*

Maximum integer values for clock settings

**THIRTY\_TWO\_BIT = 4294967295**

**class CalMicroAmpere**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *Enum*

Pre-defined microampere (uA) values for calibration current source query.

**TEN = '10uA'**

**TWO\_HUNDRED = '200uA'**

**class CjcType**(*value=<no\_arg>*, *names=None*, *module=None*, *qualname=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *NameEnum*

CJC slope and offset

```
internal = (1, 0)

lm34 = (55.56, 255.37)

class ClockFrequency(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                     start=1, boundary=None)
    Bases: IntEnum
    Available clock frequencies, in Hz
    FIVE_MHZ = 5000000
    FORTY_MHZ = 40000000
    MAXIMUM = 80000000
    MINIMUM = 312500
    TEN_MHZ = 10000000
    TWELVE_HUNDRED_FIFTY_KHZ = 1250000
    TWENTY_FIVE_HUNDRED_KHZ = 2500000
    TWENTY_MHZ = 20000000

DIOChannel
    alias of TSeriesDIOChannel

class DIOStatus(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                 boundary=None)
    Bases: IntEnum
    State of a digital I/O channel.
    HIGH = 1
    LOW = 0

class DeviceType(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                 boundary=None)
    Bases: AutoNumberNameEnum
    LabJack device types.
    Can be also looked up by ambiguous Product ID (p_id) or by instance name: `python
    LabJackDeviceType(4) is LabJackDeviceType('T4')`  

    ANY = 1
    T4 = 2
    T7 = 3
    T7_PRO = 4

    classmethod get_by_p_id(p_id: int) → Union[DeviceType, List[DeviceType]]
        Get LabJack device type instance via LabJack product ID.
        Note: Product ID is not unambiguous for LabJack devices.
```

**Parameters**

**p\_id** – Product ID of a LabJack device

**Returns**

Instance or list of instances of *LabJackDeviceType*

**Raises**

**ValueError** – when Product ID is unknown

```
class TemperatureUnit(value=<no_arg>, names=None, module=None, qualname=None, type=None,  
start=1, boundary=None)
```

Bases: [NameEnum](#)

Temperature unit (to be returned)

**C = 1**

**F = 2**

**K = 0**

```
class ThermocoupleType(value=<no_arg>, names=None, module=None, qualname=None, type=None,  
start=1, boundary=None)
```

Bases: [NameEnum](#)

Thermocouple type; NONE means disable thermocouple mode.

**C = 30**

**E = 20**

**J = 21**

**K = 22**

**NONE = 0**

**PT100 = 40**

**PT1000 = 42**

**PT500 = 41**

**R = 23**

**S = 25**

**T = 24**

```
config_high_pulse(address: Union[str, TSeriesDIOChannel], t_start: Union[int, float], t_width:  
Union[int, float], n_pulses: int = 1) → None
```

Configures one FIO channel to send a timed HIGH pulse. Configure multiple channels to send pulses with relative timing accuracy. Times have a maximum resolution of 1e-7 seconds @ 10 MHz. :param address: FIO channel: [T7] FIO0;2;3;4;5. [T4] FIO6;7. :raises LabJackError if address is not supported. :param t\_start: pulse start time in seconds. :raises ValueError: if t\_start is negative or would exceed the clock period. :param t\_width: duration of high pulse, in seconds. :raises ValueError: if t\_width is negative or would exceed the clock period. :param n\_pulses: number of pulses to be sent; single pulse default. :raises TypeError if n\_pulses is not of type int. :raises ValueError if n\_pulses is negative or exceeds the 32bit limit.

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**disable\_pulses(\*addresses: Optional[Union[str, TSeriesDIOChannel]]) → None**

Disable previously configured pulse channels. :param addresses: tuple of FIO addresses. All channels disabled if no argument is passed.

**enable\_clock(clock\_enabled: bool) → None**

Enable/disable LabJack clock to configure or send pulses. :param clock\_enabled: True -> enable, False -> disable. :raises TypeError: if clock\_enabled is not of type bool

**get\_ain(\*channels: int) → Union[float, Sequence[float]]**

Read currently measured value (voltage, resistance, ...) from one or more of analog inputs.

**Parameters**

**channels** – AIN number or numbers (0..254)

**Returns**

the read value (voltage, resistance, ...) as *float* or *tuple* of them in case multiple channels given

**get\_cal\_current\_source(name: Union[str, CalMicroAmpere]) → float**

This function will return the calibration of the chosen current source, this ist not a measurement!

The value was stored during fabrication.

**Parameters**

**name** – ‘200uA’ or ‘10uA’ current source

**Returns**

calibration of the chosen current source in ampere

**get\_clock() → Dict[str, Union[int, float]]**

Return clock settings read from LabJack.

**get\_digital\_input(address: Union[str, TSeriesDIOChannel]) → DIOStatus**

Get the value of a digital input.

allowed names for T7 (Pro): FIO0 - FIO7, EIO0 - EIO 7, CIO0- CIO3, MIO0 - MIO2 :param address: name of the output -> ‘FIO0’ :return: HIGH when *address* DIO is high, and LOW when *address* DIO is low

**get\_product\_id() → int**

This function returns the product ID reported by the connected device.

Attention: returns 7 for both T7 and T7-Pro devices!

**Returns**

integer product ID of the device

**get\_product\_name(force\_query\_id=False) → str**

This function will return the product name based on product ID reported by the device.

Attention: returns “T7” for both T7 and T7-Pro devices!

**Parameters**

**force\_query\_id** – boolean flag to force *get\_product\_id* query to device instead of using cached device type from previous queries.

**Returns**

device name string, compatible with *LabJack.DeviceType*

**get\_product\_type**(*force\_query\_id*: bool = False) → *DeviceType*

This function will return the device type based on reported device type and in case of unambiguity based on configuration of device's communication protocol (e.g. for "T7" and "T7\_PRO" devices), or, if not available first matching.

**Parameters**

**force\_query\_id** – boolean flag to force *get\_product\_id* query to device instead of using cached device type from previous queries.

**Returns**

*DeviceType* instance

**Raises**

*LabJackIdentifierDIOError* – when read Product ID is unknown

**get\_sbus\_rh**(*number*: int) → float

Read the relative humidity value from a serial SBUS sensor.

**Parameters**

**number** – port number (0..22)

**Returns**

relative humidity in %RH

**get\_sbus\_temp**(*number*: int) → float

Read the temperature value from a serial SBUS sensor.

**Parameters**

**number** – port number (0..22)

**Returns**

temperature in Kelvin

**get\_serial\_number**() → int

Returns the serial number of the connected LabJack.

**Returns**

Serial number.

**read\_resistance**(*channel*: int) → float

Read resistance from specified channel.

**Parameters**

**channel** – channel with resistor

**Returns**

resistance value with 2 decimal places

**read\_thermocouple**(*pos\_channel*: int) → float

Read the temperature of a connected thermocouple.

**Parameters**

**pos\_channel** – is the AIN number of the positive pin

**Returns**

temperature in specified unit

**send\_pulses**(\*addresses: Union[str, TSeriesDIOChannel]) → [None](#)

Sends pre-configured pulses for specified addresses. :param addresses: tuple of FIO addresses :raises LabJackError if an address has not been configured.

**set\_ain\_differential**(pos\_channel: int, differential: bool) → [None](#)

Sets an analog input to differential mode or not. T7-specific: For base differential channels, positive must be even channel from 0-12 and negative must be positive+1. For extended channels 16-127, see Mux80 datasheet.

#### Parameters

- **pos\_channel** – is the AIN number (0..12)
- **differential** – True or False

#### Raises

[LabJackError](#) – if parameters are unsupported

**set\_ain\_range**(channel: int, vrange: Union[Real, AInRange]) → [None](#)

Set the range of an analog input port.

#### Parameters

- **channel** – is the AIN number (0..254)
- **vrange** – is the voltage range to be set

**set\_ain\_resistance**(channel: int, vrange: Union[Real, AInRange], resolution: int) → [None](#)

Set the specified channel to resistance mode. It utilized the 200uA current source of the LabJack.

#### Parameters

- **channel** – channel that should measure the resistance
- **vrange** – voltage range of the channel
- **resolution** – resolution index of the channel T4: 0-5, T7: 0-8, T7-Pro 0-12

**set\_ain\_resolution**(channel: int, resolution: int) → [None](#)

Set the resolution index of an analog input port.

#### Parameters

- **channel** – is the AIN number (0..254)
- **resolution** – is the resolution index within 0...`get\_product\_type().ain\_max\_resolution` range; 0 will set the resolution index to default value.

**set\_ain\_thermocouple**(pos\_channel: int, thermocouple: Union[None, str, ThermocoupleType], cjc\_address: int = 60050, cjc\_type: Union[str, CjcType] = CjcType.internal, vrange: Union[Real, AInRange] = AInRange.ONE\_HUNDREDTH, resolution: int = 10, unit: Union[str, TemperatureUnit] = TemperatureUnit.K) → [None](#)

Set the analog input channel to thermocouple mode.

#### Parameters

- **pos\_channel** – is the analog input channel of the positive part of the differential pair
- **thermocouple** – None to disable thermocouple mode, or string specifying the thermocouple type
- **cjc\_address** – modbus register address to read the CJC temperature
- **cjc\_type** – determines cjc slope and offset, ‘internal’ or ‘lm34’

- **vrange** – measurement voltage range
- **resolution** – resolution index (T7-Pro: 0-12)
- **unit** – is the temperature unit to be returned ('K', 'C' or 'F')

**Raises**

**LabJackError** – if parameters are unsupported

**set\_analog\_output**(channel: int, value: Union[int, float]) → *None*

Set the voltage of a analog output port

**Parameters**

- **channel** – DAC channel number 1/0
- **value** – The output voltage value 0-5 Volts int/float

**set\_clock**(clock\_frequency: Union[Number, ClockFrequency] = 10000000, clock\_period: Number = 1) → *None*

Configure LabJack clock for pulse out feature. :param clock\_frequency: clock frequency in Hz; default 10 MHz for base 10. :raises ValueError: if clock\_frequency is not allowed (see ClockFrequency). :param clock\_period: clock roll time in seconds; default 1s, 0 for max. :raises ValueError: if clock\_period exceeds the 32bit tick limit. Clock period determines pulse spacing when using multi-pulse settings. Ensure period exceeds maximum intended pulse end time.

**set\_digital\_output**(address: str, state: Union[int, DIOStatus]) → *None*

Set the value of a digital output.

**Parameters**

- **address** – name of the output -> 'FIO0'
- **state** – state of the output -> *DIOStatus* instance or corresponding *int* value

**start()** → *None*

Start the Device.

**stop()** → *None*

Stop the Device.

**exception LabJackError**

Bases: *DeviceError*


General Error for the LabJack device.

**exception LabJackIdentifierDIOError**

Bases: *LabJackError*

Error indicating a wrong DIO identifier

## `hvl_ccb.dev.lauda`

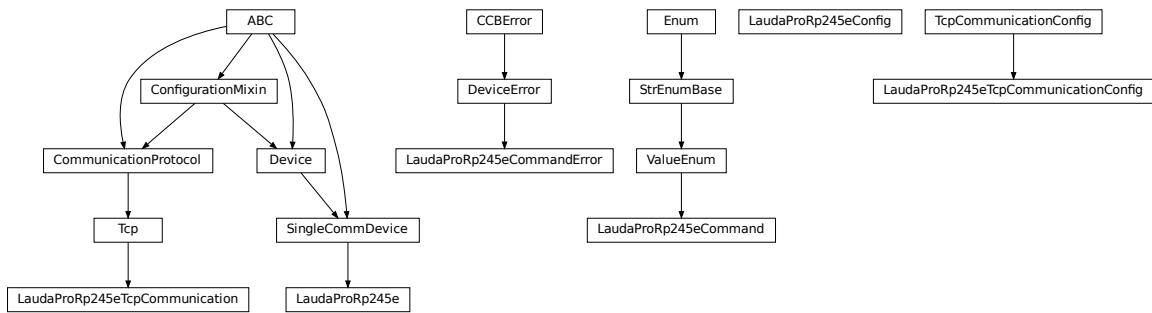

Device class for controlling a Lauda PRO RP245E, circulation chiller over TCP.

**class** `LaudaProRp245e`(*com, dev\_config=None*)

Bases: `SingleCommDevice`

Lauda RP245E circulation chiller class.

**static config\_cls()** → Type[`LaudaProRp245eConfig`]

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**continue\_ramp()** → str

Continue current ramp program.

**Returns**

reply of the device to the last call of “query”

**static default\_com\_cls()** → Type[`LaudaProRp245eTcpCommunication`]

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**get\_bath\_temp()** → float

:return : float value of measured lauda bath temp in °C

**get\_device\_type()** → str

:return : Connected Lauda device type (for connection/com test)

**pause()** → str

Stop temperature control and pump.

**Returns**

reply of the device to the last call of “query”

**pause\_ramp()** → str

Pause current ramp program.

**Returns**

reply of the device to the last call of “query”

**reset\_ramp()** → str

Delete all segments from current ramp program.

**Returns**

reply of the device to the last call of “query”

**run()** → str

Start temperature control & pump.

**Returns**

reply of the device to the last call of “query”

**set\_control\_mode(mod: Union[int, ExtControlModeEnum] = ExtControlModeEnum.INTERNAL)** → str

Define control mode. 0 = INTERNAL (control bath temp), 1 = EXPT100 (pt100 attached to chiller), 2 = ANALOG, 3 = SERIAL, 4 = USB, 5 = ETH (to be used when passing the ext. temp. via ethernet) (temperature then needs to be passed every second, when not using options 3, 4, or 5)

**Parameters**

**mod** – temp control mode (control internal temp or external temp).

**Returns**

reply of the device to the last call of “query” (“OK”, if command was recognized”)

**set\_external\_temp(external\_temp: float = 20.0)** → str

Pass value of external controlled temperature. Should be done every second, when control of external temperature is active. Has to be done right before control of external temperature is activated.

**Parameters**

**external\_temp** – current value of external temperature to be controlled.

**Returns**

reply of the device to the last call of “query”

**set\_pump\_level(pump\_level: int = 6)** → str

Set pump level Raises ValueError, if pump level is invalid.

**Parameters**

**pump\_level** – pump level.

**Returns**

reply of the device to the last call of “query”

**set\_ramp\_iterations(num: int = 1)** → str

Define number of ramp program cycles.

**Parameters**

**num** – number of program cycles to be performed.

**Returns**

reply of the device to the last call of “query”

**set\_ramp\_program(program: int = 1)** → str

Define ramp program for following ramp commands. Raises ValueError if maximum number of ramp programs (5) is exceeded.

**Parameters**

**program** – Number of ramp program to be activated for following commands.

**Returns**

reply of the device to the last call of “query”

**set\_ramp\_segment**(*temp: float = 20.0, dur: int = 0, tol: float = 0.0, pump: int = 6*) → str

Define segment of current ramp program - will be attached to current program. Raises ValueError, if pump level is invalid.

**Parameters**

- **temp** – target temperature of current ramp segment
- **dur** – duration in minutes, in which target temperature should be reached
- **tol** – tolerance at which target temperature should be reached (for 0.00, next segment is started after dur has passed).
- **pump** – pump level to be used for this program segment.

**Returns**

reply of the device to the last call of “query”

**set\_temp\_set\_point**(*temp\_set\_point: float = 20.0*) → str

Define temperature set point

**Parameters**

**temp\_set\_point** – temperature set point.

**Returns**

reply of the device to the last call of “query”

**start()** → *None*

Start this device.

**start\_ramp()** → str

Start current ramp program.

**Returns**

reply of the device to the last call of “query”

**stop()** → *None*

Stop this device. Disables access and closes the communication protocol.

**stop\_ramp()** → str

Stop current ramp program.

**Returns**

reply of the device to the last call of “query”

**validate\_pump\_level**(*level: int*)

Validates pump level. Raises ValueError, if pump level is incorrect. :param level: pump level, integer

**class LaudaProRp245eCommand**(*value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None*)

Bases: *ValueEnum*

Commands for Lauda PRO RP245E Chiller Command strings most often need to be complimented with a parameter (attached as a string) before being sent to the device. Commands implemented as defined in “Lauda Betriebsanleitung fuer PRO Badthermostate und Umwaelzthermostate” pages 42 - 49

**BATH\_TEMP = 'IN\_PV\_00'**

Request internal bath temperature

**COM\_TIME\_OUT = 'OUT\_SP\_08\_'**

Define communication time out

---

```

CONT_MODE = 'OUT_MODE_01_'
    Set control mode 1=internal, 2=ext. analog, 3=ext. serial, 4=USB, 5=ethernet

DEVICE_TYPE = 'TYPE'
    Request device type

EXTERNAL_TEMP = 'OUT_PV_05_'
    Pass on external controlled temperature

LOWER_TEMP = 'OUT_SP_05_'
    Define lower temp limit

OPERATION_MODE = 'OUT_SP_02_'
    Define operation mode

PUMP_LEVEL = 'OUT_SP_01_'
    Define pump level 1-8

RAMP_CONTINUE = 'RMP_CONT'
    Continue a paused ramp program

RAMP_DELETE = 'RMP_RESET'
    Reset a selected ramp program

RAMP_ITERATIONS = 'RMP_OUT_02_'
    Define how often a ramp program should be iterated

RAMP_PAUSE = 'RMP_PAUSE'
    Pause a selected ramp program

RAMP_SELECT = 'RMP_SELECT_'
    Select a ramp program (target for all further ramp commands)

RAMP_SET = 'RMP_OUT_00_'
    Define parameters of a selected ramp program

RAMP_START = 'RMP_START'
    Start a selected ramp program

RAMP_STOP = 'RMP_STOP'
    Stop a running ramp program

START = 'START'
    Start temp control (pump and heating/cooling)

STOP = 'STOP'
    Stop temp control (pump and heating/cooling)

TEMP_SET_POINT = 'OUT_SP_00_'
    Define temperature set point

UPPER_TEMP = 'OUT_SP_04_'
    Define upper temp limit

build_str(param: str = "", terminator: str = '\r\n')
    Build a command string for sending to the device

```

#### Parameters

- **param** – Command's parameter given as string

- **terminator** – Command's terminator

**Returns**

Command's string with a parameter and terminator

**exception LaudaProRp245eCommandError**

Bases: *DeviceError*

Error raised when an error is returned upon a command.

```
class LaudaProRp245eConfig(temp_set_point_init: Union[int, float] = 20.0, pump_init: int = 6, upper_temp:  
    Union[int, float] = 80.0, lower_temp: Union[int, float] = -55.0, com_time_out:  
    Union[int, float] = 0, max_pump_level: int = 8, max_pr_number: int = 5,  
    operation_mode: Union[int, OperationModeEnum] =  
        OperationModeEnum.AUTO, control_mode: Union[int, ExtControlModeEnum]  
        = ExtControlModeEnum.INTERNAL)
```

Bases: *object*

Configuration for the Lauda RP245E circulation chiller.

**class ExtControlModeEnum(*value*)**

Bases: *IntEnum*

Source for definition of external, controlled temperature (option 2, 3 and 4 are not available with current configuration of the Lauda RP245E, add-on hardware would required)

**ANALOG = 2**

**ETH = 5**

**EXPT100 = 1**

**INTERNAL = 0**

**SERIAL = 3**

**USB = 4**

**class OperationModeEnum(*value*)**

Bases: *IntEnum*

Operation Mode (Cooling OFF/Cooling On/AUTO - set to AUTO)

**AUTO = 2**

Automatically select heating/cooling

**COOLOFF = 0**

**COOLON = 1**

**clean\_values() → *None***

**com\_time\_out: Union[int, float] = 0**

Communication time out (0 = OFF)

**control\_mode: Union[int, ExtControlModeEnum] = 0**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True****classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**lower\_temp: Union[int, float] = -55.0**

Lower temperature limit (safe for Galden HT135 cooling liquid)

**max\_pr\_number: int = 5**

Maximum number of ramp programs that can be stored in the memory of the chiller

**max\_pump\_level: int = 8**

Highest pump level of the chiller

**operation\_mode: Union[int, OperationModeEnum] = 2****classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**pump\_init: int = 6**

Default pump Level

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**temp\_set\_point\_init: Union[int, float] = 20.0**

Default temperature set point

**upper\_temp: Union[int, float] = 80.0**

Upper temperature limit (safe for Galden HT135 cooling liquid)

**class LaudaProRp245eTcpCommunication(*configuration*)**

Bases: [Tcp](#)

Implements the Communication Protocol for Lauda PRO RP245E TCP connection.

**close()** → *None*

Close the Lauda PRO RP245E TCP connection.

**static config\_cls()** → Type[*LaudaProRp245eTcpCommunicationConfig*]

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**open()** → *None*

Open the Lauda PRO RP245E TCP connection.

**Raises**

*LaudaProRp245eCommandError* – if the connection fails.

**query\_command(command: LaudaProRp245eCommand, param: str = "")** → str

Send and receive function. E.g. to be used when setting/changing device setting. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None

**read()** → str

Receive value function. :return: reply from device as a string, the terminator, as well as the 'OK' stripped from the reply to make it directly useful as a value (e.g. in case the internal bath temperature is requested)

**write\_command(command: LaudaProRp245eCommand, param: str = "")** → *None*

Send command function. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None

**class LaudaProRp245eTcpCommunicationConfig(host: Union[str, IPv4Address, IPv6Address], port: int = 54321, bufsize: int = 1024, wait\_sec\_pre\_read\_or\_write: Union[int, float] = 0.005, terminator: str = '\r\n')**

Bases: *TcpCommunicationConfig*

Configuration dataclass for *LaudaProRp245eTcpCommunication*.

**clean\_values()** → *None*

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.


```
terminator: str = '\r\n'
```

The terminator character

```
wait_sec_pre_read_or_write: Union[int, float] = 0.005
```

Delay time between commands in seconds

## hvl\_ccb.dev.mbw973

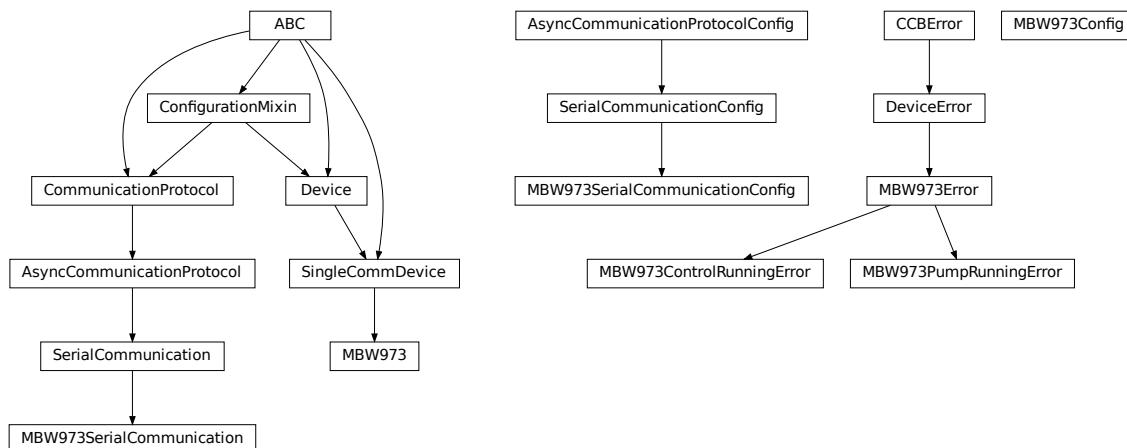

Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

```
class MBW973(com, dev_config=None)
```

Bases: *SingleCommDevice*

MBW 973 dew point mirror device class.

```
static config_cls()
```

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**is\_done()** → bool

Poll status of the dew point mirror and return True, if all measurements are done.

**Returns**

True, if all measurements are done; False otherwise.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**read(*cast\_type*: ~typing.Type = <class 'str'>)**

Read value from *self.com* and cast to *cast\_type*. Raises *ValueError* if read text (*str*) is not convertible to *cast\_type*, e.g. to *float* or to *int*.

**Returns**

Read value of *cast\_type* type.

**read\_float()** → float

Convenience wrapper for *self.read()*, with typing hint for return value.

**Returns**

Read *float* value.

**read\_int()** → int

Convenience wrapper for *self.read()*, with typing hint for return value.

**Returns**

Read *int* value.

**read\_measurements()** → Dict[str, float]

Read out measurement values and return them as a dictionary.

**Returns**

Dictionary with values.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**set\_measuring\_options(*humidity*: bool = True, *sf6\_purity*: bool = False) → None**

Send measuring options to the dew point mirror.

**Parameters**

- **humidity** – Perform humidity test or not?
- **sf6\_purity** – Perform SF6 purity test or not?

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**start()** → *None*

Start this device. Opens the communication protocol and retrieves the set measurement options from the device.

**Raises**

*SerialCommunicationIOError* – when communication port cannot be opened.

**start\_control()** → *None*

Start dew point control to acquire a new value set.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**stop()** → *None*

Stop the device. Closes also the communication protocol.

**write(*value*)** → *None*

Send *value* to *self.com*.

**Parameters**

**value** – Value to send, converted to *str*.

**Raises**

*SerialCommunicationIOError* – when communication port is not opened

**class MBW973Config(*polling\_interval*: Union[int, float] = 2)**

Bases: *object*

Device configuration dataclass for MBW973.

**clean\_values()****force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True****classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**polling\_interval: Union[int, float] = 2**

Polling period for *is\_done* status queries [in seconds].

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**exception MBW973ControlRunningError**

Bases: *MBW973Error*

Error indicating there is still a measurement running, and a new one cannot be started.

**exception MBW973Error**

Bases: *DeviceError*

General error with the MBW973 dew point mirror device.

**exception MBW973PumpRunningError**

Bases: *MBW973Error*

Error indicating the pump of the dew point mirror is still recovering gas, unable to start a new measurement.

**class MBW973SerialCommunication(*configuration*)**

Bases: *SerialCommunication*

Specific communication protocol implementation for the MBW973 dew point mirror. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class MBW973SerialCommunicationConfig(*terminator: bytes = b'\r', encoding: str = 'utf-8', encoding\_error\_handling: str = 'strict', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3*)**

Bases: *SerialCommunicationConfig*

**baudrate: int = 9600**

Baudrate for MBW973 is 9600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

MBW973 does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**stopbits: Union[int, SerialCommunicationStopbits] = 1**

MBW973 does use one stop bit


**terminator: bytes = b'\r'**

The terminator is only CR

**timeout: Union[int, float] = 3**

use 3 seconds timeout as default

## hvl\_ccb.dev.newport

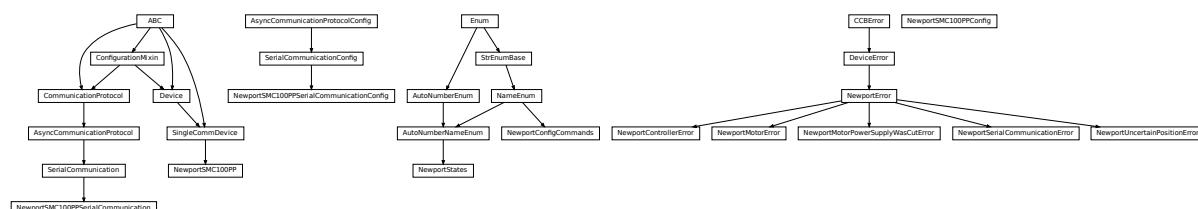

Device class for Newport SMC100PP stepper motor controller with serial communication.

The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/f/smci00-single-axis-dc-or-stepper-motion-controller>

**class NewportConfigCommands(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *NameEnum*

Commands predefined by the communication protocol of the SMC100PP

**AC = 'acceleration'**

**BA = 'backlash\_compensation'**

**BH = 'hysteresis\_compensation'**

**FRM = 'micro\_step\_per\_full\_step\_factor'**

```
FRS = 'motion_distance_per_full_step'
HT = 'home_search_type'
JR = 'jerk_time'
OH = 'home_search_velocity'
OT = 'home_search_timeout'
QIL = 'peak_output_current_limit'
SA = 'rs485_address'
SL = 'negative_software_limit'
SR = 'positive_software_limit'
VA = 'velocity'
VB = 'base_velocity'
ZX = 'stage_configuration'

exception NewportControllerError
    Bases: NewportError
    Error with the Newport controller.

exception NewportError
    Bases: DeviceError
    General Error for Newport Device

exception NewportMotorError
    Bases: NewportError
    Error with the Newport motor.

exception NewportMotorPowerSupplyWasCutError
    Bases: NewportError
    Error with the Newport motor after the power supply was cut and then restored, without interrupting the communication with the controller.

class NewportSMC100PP(com, dev_config=None)
    Bases: SingleCommDevice
    Device class of the Newport motor controller SMC100PP

    class MotorErrors(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
        Bases: Enum
        Possible motor errors reported by the motor during get_state().
        DC_VOLTAGE_TOO_LOW = 3
        FOLLOWING_ERROR = 6
        HOMING_TIMEOUT = 5
```

```
NED_END_OF_TURN = 11
OUTPUT_POWER_EXCEEDED = 2
PEAK_CURRENT_LIMIT = 9
POS_END_OF_TURN = 10
RMS_CURRENT_LIMIT = 8
SHORT_CIRCUIT = 7
WRONG_ESP_STAGE = 4

class StateMessages(value=<no_arg>, names=None, module=None, qualname=None, type=None,
                    start=1, boundary=None)
Bases: Enum
Possible messages returned by the controller on get_state() query.

CONFIG = '14'

DISABLE_FROM_JOGGING = '3E'
DISABLE_FROM_MOVING = '3D'
DISABLE_FROM_READY = '3C'
HOMING_FROM_RS232 = '1E'
HOMING_FROM_SMC = '1F'
JOGGING_FROM_DISABLE = '47'
JOGGING_FROM_READY = '46'
MOVING = '28'
NO_REF_ESP_STAGE_ERROR = '10'
NO_REF_FROM_CONFIG = '0C'
NO_REF_FROM_DISABLED = '0D'
NO_REF_FROM_HOMING = '0B'
NO_REF_FROM_JOGGING = '11'
NO_REF_FROM_MOVING = '0F'
NO_REF_FROM_READY = '0E'
NO_REF_FROM_RESET = '0A'
READY_FROM_DISABLE = '34'
READY_FROM_HOMING = '32'
READY_FROM_JOGGING = '35'
READY_FROM_MOVING = '33'
```

**States**

alias of [NewportStates](#)

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**exit\_configuration(*add: Optional[int] = None*) → *None***

Exit the CONFIGURATION state and go back to the NOT REFERENCED state. All configuration parameters are saved to the device's memory.

**Parameters**

**add** – controller address (1 to 31)

**Raises**

- [SerialCommunicationIOError](#) – if the com is closed
- [NewportSerialCommunicationError](#) – if an unexpected answer is obtained
- [NewportControllerError](#) – if the controller reports an error

**get\_acceleration(*add: Optional[int] = None*) → Union[int, float]**

Leave the configuration state. The configuration parameters are saved to the device's memory.

**Parameters**

**add** – controller address (1 to 31)

**Returns**

acceleration (preset units/s<sup>2</sup>), value between 1e-6 and 1e12

**Raises**

- [SerialCommunicationIOError](#) – if the com is closed
- [NewportSerialCommunicationError](#) – if an unexpected answer is obtained
- [NewportControllerError](#) – if the controller reports an error

**get\_controller\_information(*add: Optional[int] = None*) → str**

Get information on the controller name and driver version

**Parameters**

**add** – controller address (1 to 31)

**Returns**

controller information

**Raises**

- [SerialCommunicationIOError](#) – if the com is closed
- [NewportSerialCommunicationError](#) – if an unexpected answer is obtained
- [NewportControllerError](#) – if the controller reports an error

**get\_motor\_configuration**(*add: Optional[int] = None*) → Dict[str, float]

Query the motor configuration and returns it in a dictionary.

**Parameters**

**add** – controller address (1 to 31)

**Returns**

dictionary containing the motor's configuration

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**get\_move\_duration**(*dist: Union[int, float], add: Optional[int] = None*) → float

Estimate the time necessary to move the motor of the specified distance.

**Parameters**

- **dist** – distance to travel
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**get\_negative\_software\_limit**(*add: Optional[int] = None*) → Union[int, float]

Get the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

**Parameters**

**add** – controller address (1 to 31)

**Returns**

negative software limit (preset units), value between -1e12 and 0

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**get\_position**(*add: Optional[int] = None*) → float

Returns the value of the current position.

**Parameters**

**add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error
- **NewportUncertainPositionError** – if the position is ambiguous

**get\_positive\_software\_limit**(*add: Optional[int] = None*) → Union[int, float]

Get the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

**Parameters**

**add** – controller address (1 to 31)

**Returns**

positive software limit (preset units), value between 0 and 1e12

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**get\_state**(*add: int = None*) → *StateMessages*

Check on the motor errors and the controller state

**Parameters**

**add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error
- **NewportMotorError** – if the motor reports an error

**Returns**

state message from the device (member of StateMessages)

**go\_home**(*add: Optional[int] = None*) → *None*

Move the motor to its home position.

**Parameters**

**add** – controller address (1 to 31), defaults to self.address

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**go\_to\_configuration**(*add: Optional[int] = None*) → *None*

This method is executed during start(). It can also be executed after a reset(). The controller is put in CONFIG state, where configuration parameters can be changed.

**Parameters**

**add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**initialize**(*add: Optional[int] = None*) → *None*

Puts the controller from the NOT\_REF state to the READY state. Sends the motor to its “home” position.

**Parameters**

**add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**move\_to\_absolute\_position**(*pos: Union[int, float], add: Optional[int] = None*) → *None*

Move the motor to the specified position.

**Parameters**

- **pos** – target absolute position (affected by the configured offset)
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**move\_to\_relative\_position**(*pos: Union[int, float], add: Optional[int] = None*) → *None*

Move the motor of the specified distance.

**Parameters**

- **pos** – distance to travel (the sign gives the direction)
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**reset**(*add: Optional[int] = None*) → *None*

Resets the controller, equivalent to a power-up. This puts the controller back to NOT REFERENCED state, which is necessary for configuring the controller.

**Parameters**

**add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**set\_acceleration**(*acc: Union[int, float], add: Optional[int] = None*) → *None*

Leave the configuration state. The configuration parameters are saved to the device’s memory.

**Parameters**

- **acc** – acceleration (preset units/s<sup>2</sup>), value between 1e-6 and 1e12
- **add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**set\_motor\_configuration**(*add*: *Optional[int]* = *None*, *config*: *Optional[dict]* = *None*) → *None*

Set the motor configuration. The motor must be in CONFIG state.

**Parameters**

- **add** – controller address (1 to 31)
- **config** – dictionary containing the motor's configuration

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**set\_negative\_software\_limit**(*lim*: *Union[int, float]*, *add*: *Optional[int]* = *None*) → *None*

Set the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

**Parameters**

- **lim** – negative software limit (preset units), value between -1e12 and 0
- **add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**set\_positive\_software\_limit**(*lim*: *Union[int, float]*, *add*: *Optional[int]* = *None*) → *None*

Set the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

**Parameters**

- **lim** – positive software limit (preset units), value between 0 and 1e12
- **add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**start()**

Opens the communication protocol and applies the config.

**Raises**

- `SerialCommunicationIOError` – when communication port cannot be opened

**stop() → None**

Stop the device. Close the communication protocol.

**stop\_motion(add: Optional[int] = None) → None**

Stop a move in progress by decelerating the positioner immediately with the configured acceleration until it stops. If a controller address is provided, stops a move in progress on this controller, else stops the moves on all controllers.

**Parameters**

`add` – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**wait\_until\_motor\_initialized(add: Optional[int] = None) → None**

Wait until the motor leaves the HOMING state (at which point it should have arrived to the home position).

**Parameters**

`add` – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

```
class NewportSMC100PPConfig(address: int = 1, user_position_offset: Union[int, float] = 23.987,
                             screw_scaling: Union[int, float] = 1, exit_configuration_wait_sec: Union[int, float] = 5, move_wait_sec: Union[int, float] = 1, acceleration: Union[int, float] = 10, backlash_compensation: Union[int, float] = 0, hysteresis_compensation: Union[int, float] = 0.015, micro_step_per_full_step_factor: int = 100, motion_distance_per_full_step: Union[int, float] = 0.01, home_search_type: Union[int, HomeSearch] = HomeSearch.HomeSwitch, jerk_time: Union[int, float] = 0.04, home_search_velocity: Union[int, float] = 4, home_search_timeout: Union[int, float] = 27.5, home_search_polling_interval: Union[int, float] = 1, peak_output_current_limit: Union[int, float] = 0.4, rs485_address: int = 2, negative_software_limit: Union[int, float] = -23.5, positive_software_limit: Union[int, float] = 25, velocity: Union[int, float] = 4, base_velocity: Union[int, float] = 0, stage_configuration: Union[int, EspStageConfig] = EspStageConfig.EnableEspStageCheck)
```

Bases: `object`

Configuration dataclass for the Newport motor controller SMC100PP.

```
class EspStageConfig(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: `IntEnum`

Different configurations to check or not the motor configuration upon power-up.

```
DisableEspStageCheck = 1
EnableEspStageCheck = 3
UpdateEspStageInfo = 2

class HomeSearch(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
Bases: IntEnum
Different methods for the motor to search its home position during initialization.

CurrentPosition = 1
EndOfRunSwitch = 4
EndOfRunSwitch_and_Index = 3
HomeSwitch = 2
HomeSwitch_and_Index = 0

acceleration: Union[int, float] = 10
address: int = 1
backlash_compensation: Union[int, float] = 0
base_velocity: Union[int, float] = 0
clean_values()
exit_configuration_wait_sec: Union[int, float] = 5
force_value(fieldname, value)
Forces a value to a dataclass field despite the class being frozen.
NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.
```

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

```
home_search_polling_interval: Union[int, float] = 1
home_search_timeout: Union[int, float] = 27.5
home_search_type: Union[int, HomeSearch] = 2
home_search_velocity: Union[int, float] = 4
hysteresis_compensation: Union[int, float] = 0.015
is_configdataclass = True
jerk_time: Union[int, float] = 0.04
```

```
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns
        a list of strings containing all keys.

micro_step_per_full_step_factor: int = 100
motion_distance_per_full_step: Union[int, float] = 0.01
property motor_config: Dict[str, float]
    Gather the configuration parameters of the motor into a dictionary.

    Returns
        dict containing the configuration parameters of the motor

move_wait_sec: Union[int, float] = 1
negative_software_limit: Union[int, float] = -23.5
classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns
        a list of strings containing all optional keys.

peak_output_current_limit: Union[int, float] = 0.4
positive_software_limit: Union[int, float] = 25
post_force_value(fieldname, value)

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns
        a list of strings containing all required keys.

rs485_address: int = 2
screw_scaling: Union[int, float] = 1
stage_configuration: Union[int, EspStageConfig] = 3
user_position_offset: Union[int, float] = 23.987
velocity: Union[int, float] = 4

class NewportSMC100PPSSerialCommunication(configuration)
    Bases: SerialCommunication
    Specific communication protocol implementation for NewportSMC100 controller. Already predefines device-specific protocol parameters in config.

class ControllerErrors(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
    Bases: Enum
    Possible controller errors with values as returned by the device in response to sent commands.
```

```
ADDR_INCORRECT = 'B'  
CMD_EXEC_ERROR = 'V'  
CMD_NOT_ALLOWED = 'D'  
CMD_NOT_ALLOWED_CC = 'X'  
CMD_NOT_ALLOWED_CONFIGURATION = 'I'  
CMD_NOT_ALLOWED_DISABLE = 'J'  
CMD_NOT_ALLOWED_HOMING = 'L'  
CMD_NOT_ALLOWED_MOVING = 'M'  
CMD_NOT_ALLOWED_NOT_REFERENCED = 'H'  
CMD_NOT_ALLOWED_PP = 'W'  
CMD_NOT_ALLOWED_READY = 'K'  
CODE_OR_ADDR_INVALID = 'A'  
COM_TIMEOUT = 'S'  
DISPLACEMENT_OUT_OF_LIMIT = 'G'  
EEPROM_ACCESS_ERROR = 'U'  
ESP_STAGE_NAME_INVALID = 'F'  
HOME_STARTED = 'E'  
NO_ERROR = '@'  
PARAM_MISSING_OR_INVALID = 'C'  
POSITION_OUT_OF_LIMIT = 'N'
```

**check\_for\_error(*add*: int) → *None***

Ask the Newport controller for the last error it recorded.

This method is called after every command or query.

**Parameters**

*add* – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**query**(*add*: int, *cmd*: str, *param*: Optional[Union[int, float, str]] = None) → str

Send a query to the controller, read the answer, and check for errors. The prefix *add+cmd* is removed from the answer.

#### Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

#### Returns

the answer from the device without the prefix

#### Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**query\_multiple**(*add*: int, *cmd*: str, *prefixes*: List[str]) → List[str]

Send a query to the controller, read the answers, and check for errors. The prefixes are removed from the answers.

#### Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **prefixes** – prefixes of each line expected in the answer

#### Returns

list of answers from the device without prefix

#### Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**read\_text**() → str

Read one line of text from the serial port, and check for presence of a null char which indicates that the motor power supply was cut and then restored. The input buffer may hold additional data afterwards, since only one line is read.

This method uses *self.access\_lock* to ensure thread-safety.

#### Returns

String read from the serial port; ‘’ if there was nothing to read.

#### Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **NewportMotorPowerSupplyWasCutError** – if a null char is read

**send\_command**(*add*: int, *cmd*: str, *param*: Optional[Union[int, float, str]] = None) → None

Send a command to the controller, and check for errors.

#### Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

#### Raises

- ***SerialCommunicationIOError*** – if the com is closed
- ***NewportSerialCommunicationError*** – if an unexpected answer is obtained
- ***NewportControllerError*** – if the controller reports an error

**send\_stop**(*add*: int) → *None*

Send the general stop ST command to the controller, and check for errors.

#### Parameters

**add** – the controller address (1 to 31)

#### Returns

ControllerErrors reported by Newport Controller

#### Raises

- ***SerialCommunicationIOError*** – if the com is closed
- ***NewportSerialCommunicationError*** – if an unexpected answer is obtained

```
class NewportSMC100PPSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'ascii',
                                                encoding_error_handling: str = 'replace',
                                                wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                                default_n_attempts_read_text_nonempty: int = 10,
                                                port: Union[str, NoneType] = None, baudrate: int =
57600, parity: Union[str,
hvl_ccb.comm.serial.SerialCommunicationParity] =
<SerialCommunicationParity.NONE: 'N'>, stopbits:
Union[int,
hvl_ccb.comm.serial.SerialCommunicationStopbits] =
<SerialCommunicationStopbits.ONE: 1>, bytesize:
Union[int,
hvl_ccb.comm.serial.SerialCommunicationBytesize] =
<SerialCommunicationBytesize.EIGHTBITS: 8>,
timeout: Union[int, float] = 10)
```

Bases: *SerialCommunicationConfig*

**baudrate: int = 57600**

Baudrate for NewportSMC100 controller is 57600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

NewportSMC100 controller uses 8 bits for one data byte

**encoding: str = 'ascii'**

use ASCII as de-/encoding, cf. the manual

**encoding\_error\_handling: str = 'replace'**

replace bytes with instead of raising utf-8 exception when decoding fails

**force\_value(*fieldname*, *value*)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

#### Returns

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### Returns

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

NewportSMC100 controller does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

#### Returns

a list of strings containing all required keys.

**stopbits: Union[int, SerialCommunicationStopbits] = 1**

NewportSMC100 controller uses one stop bit

**terminator: bytes = b'\r\n'**

The terminator is CR/LF

**timeout: Union[int, float] = 10**

use 10 seconds timeout as default

**exception NewportSerialCommunicationError**

Bases: `NewportError`

Communication error with the Newport controller.

**class NewportStates(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: `AutoNumberNameEnum`

States of the Newport controller. Certain commands are allowed only in certain states.

**CONFIG = 3**

**DISABLE = 6**

**HOMING = 2**

**JOGGING = 7**

**MOVING = 5**

**NO\_REF** = 1


**READY** = 4

**exception NewportUncertainPositionError**

Bases: *NewportError*

Error with the position of the Newport motor.

## **hvl\_ccb.dev.pfeiffer\_tpg**

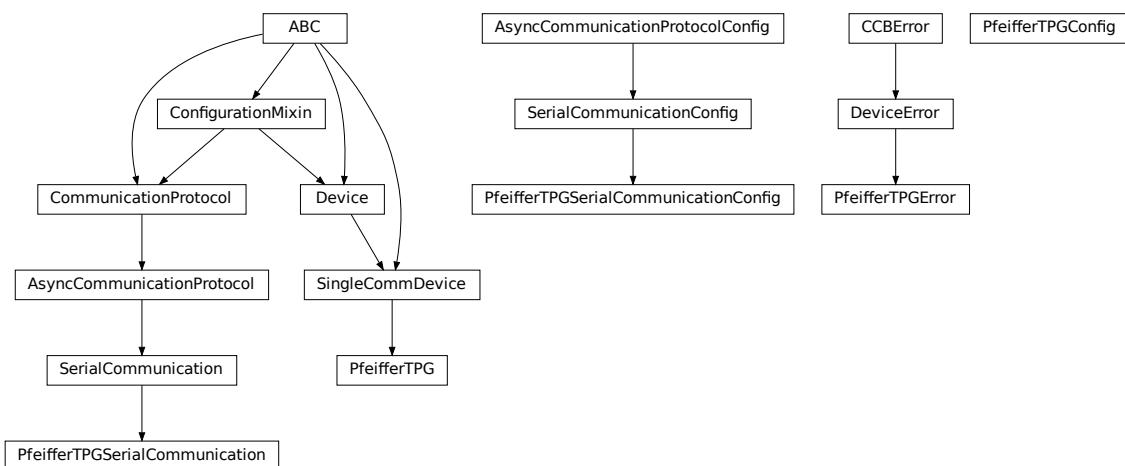

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/> measurement/activeline/controllers/

**class PfeifferTPG**(*com, dev\_config=None*)

Bases: *SingleCommDevice*

Pfeiffer TPG control unit device class

**class SensorStatus**(*value*)

Bases: *IntEnum*

An enumeration.

**Identification\_error** = 6

**No\_sensor** = 5

**Ok** = 0

**OVERRANGE** = 2

**Sensor\_error** = 3

```
Sensor_off = 4  
Underrange = 1  
class SensorTypes(value)  
    Bases: Enum  
    An enumeration.  
    CMR = 4  
    IKR = 2  
    IKR11 = 2  
    IKR9 = 2  
    IMR = 5  
    None = 7  
    PBR = 6  
    PKR = 3  
    TPR = 1  
    noSENSOR = 7  
    noSen = 7
```

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns**

the type of the standard communication protocol for this device

**get\_full\_scale\_mbar() → List[Union[int, float]]**

Get the full scale range of the attached sensors

**Returns**

full scale range values in mbar, like [0.01, 1, 0.1, 1000, 50000, 10]

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**get\_full\_scale\_unitless() → List[int]**

Get the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

**Returns**

list of full scale range values, like [0, 1, 3, 3, 2, 0]

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **PfeifferTPGError** – if command fails

**identify\_sensors()** → *None*

Send identification request TID to sensors on all channels.

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **PfeifferTPGError** – if command fails

**measure(channel: int)** → Tuple[str, float]

Get the status and measurement of one sensor

**Parameters**

**channel** – int channel on which the sensor is connected, with  $1 \leq \text{channel} \leq \text{number\_of\_sensors}$

**Returns**

measured value as float if measurement successful, sensor status as string if not

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **PfeifferTPGError** – if command fails

**measure\_all()** → List[Tuple[str, float]]

Get the status and measurement of all sensors (this command is not available on all models)

**Returns**

list of measured values as float if measurements successful, and or sensor status as strings if not

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **PfeifferTPGError** – if command fails

**property number\_of\_sensors**

**set\_full\_scale\_mbar(fsr: List[Union[int, float]])** → *None*

Set the full scale range of the attached sensors (in unit mbar)

**Parameters**

**fsr** – full scale range values in mbar, for example  $[0.01, 1000]$

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened
- **PfeifferTPGError** – if command fails

**set\_full\_scale\_unitless(fsr: List[int])** → *None*

Set the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

**Parameters**

**fsr** – list of full scale range values, like  $[0, 1, 3, 3, 2, 0]$

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened

- **PfeifferTPGError** – if command fails

**start()** → *None*

Start this device. Opens the communication protocol, and identify the sensors.

**Raises**

**SerialCommunicationIOError** – when communication port cannot be opened

**stop()** → *None*

Stop the device. Closes also the communication protocol.

**property unit**

The pressure unit of readings is always mbar, regardless of the display unit.

**class PfeifferTPGConfig(model: Union[str, Model] = Model.TPG25xA)**

Bases: object

Device configuration dataclass for Pfeiffer TPG controllers.

**class Model(value=<no\_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)**

Bases: *NameEnum*

An enumeration.

TPG25xA = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

TPGx6x = {0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8, 50000: 9}

**is\_valid\_scale\_range\_reversed\_str(v: str) → bool**

Check if given string represents a valid reversed scale range of a model.

**Parameters**

v – Reversed scale range string.

**Returns**

*True* if valid, *False* otherwise.

**clean\_values()**

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

```
model: Union[str, Model] = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4,
5000: 5, 10000: 6, 50000: 7}
```

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**exception PfeifferTPGError**

Bases: *DeviceError*

Error with the Pfeiffer TPG Controller.

**class PfeifferTPGSerialCommunication(configuration)**

Bases: *SerialCommunication*

Specific communication protocol implementation for Pfeiffer TPG controllers. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**query(cmd: str) → str**

Send a query, then read and returns the first line from the com port.

**Parameters**

**cmd** – query message to send to the device

**Returns**

first line read on the com

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if the device does not acknowledge the command or if the answer from the device is empty

**send\_command(cmd: str) → None**

Send a command to the device and check for acknowledgement.

**Parameters**

**cmd** – command to send to the device

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if the answer from the device differs from the expected acknowledgement character ‘chr(6)’.

```
class PfeifferTPGSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                             Union[str,
                                               hvl_ccb.comm.serial.SerialCommunicationParity] =
                                               <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                               Union[int,
                                                 hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                 <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                 Union[int,
                                                   hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                   <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                                   Union[int, float] = 3)
```

Bases: *SerialCommunicationConfig*

**baudrate: int = 9600**

Baudrate for Pfeiffer TPG controllers is 9600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

Pfeiffer TPG controllers do not use parity

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

```
stopbits: Union[int, SerialCommunicationStopbits] = 1
    Pfeiffer TPG controllers use one stop bit
terminator: bytes = b'\r\n'
    The terminator is <CR><LF>
timeout: Union[int, float] = 3
    use 3 seconds timeout as default
```

## hvl\_ccb.dev.picotech\_pt104


**NOTE:** PicoSDK Python wrappers already on import attempt to load the PicoSDK library; thus, the API docs can only be generated in a system with the latter installed and are by default disabled.

To build the API docs for this submodule locally edit the `docs/hvl_ccb.dev.picotech_pt104.rst` file to remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.picotech_pt104
:parts: 1

.. automodule:: hvl_ccb.dev.picotech_pt104
:members:
:undoc-members:
:show-inheritance:
```

## hvl\_ccb.dev.rs\_rto1024

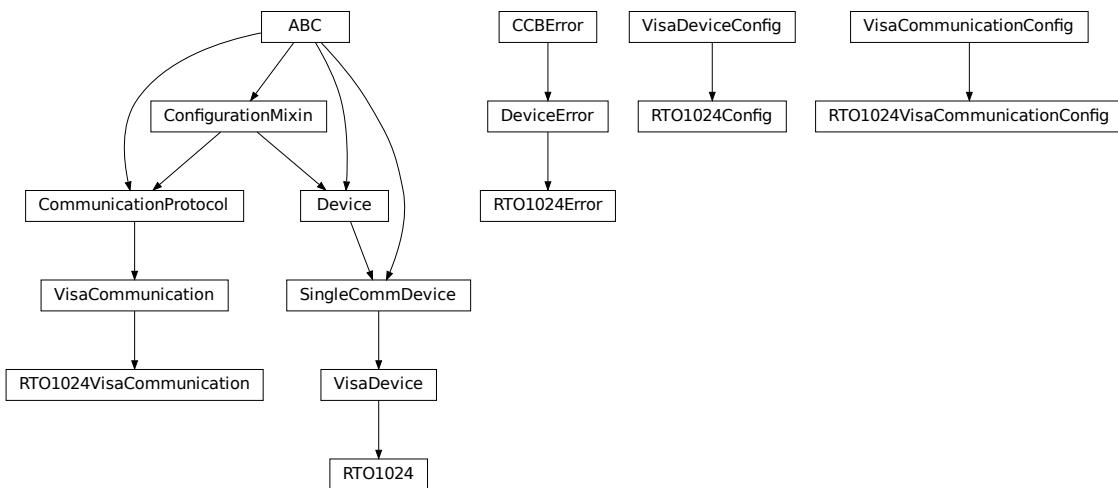

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

```
class RT01024(com: Union[RTO1024VisaCommunication, RTO1024VisaCommunicationConfig, dict],
              dev_config: Union[RTO1024Config, dict])
```

Bases: *VisaDevice*

Device class for the Rhode & Schwarz RTO 1024 oscilloscope.

**class TriggerModes**(*value*=<no\_arg>, *names*=None, *module*=None, *qualname*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *AutoNumberNameEnum*

Enumeration for the three available trigger modes.

**AUTO** = 1

**FREERUN** = 3

**NORMAL** = 2

**classmethod names()**

Returns a list of the available trigger modes. :return: list of strings

**activate\_measurements**(*meas\_n*: int, *source*: str, *measurements*: List[str], *category*: str = 'AMPTime')

Activate the list of ‘measurements’ of the waveform ‘source’ in the measurement box number ‘meas\_n’. The list ‘measurements’ starts with the main measurement and continues with additional measurements of the same ‘category’.

#### Parameters

- **meas\_n** – measurement number 1..8
- **source** – measurement source, for example C1W1
- **measurements** – list of measurements, the first one will be the main measurement.
- **category** – the category of measurements, by default AMPTime

**backup\_waveform**(*filename*: str) → *None*

Backup a waveform file from the standard directory specified in the device configuration to the standard backup destination specified in the device configuration. The filename has to be specified without .bin or path.

#### Parameters

**filename** – The waveform filename without extension and path

**static config\_cls()**

Return the default configdataclass class.

#### Returns

a reference to the default configdataclass class

**static default\_com\_cls()**

Return the default communication protocol for this device type, which is VisaCommunication.

#### Returns

the VisaCommunication class

**file\_copy**(*source*: str, *destination*: str) → *None*

Copy a file from one destination to another on the oscilloscope drive. If the destination file already exists, it is overwritten without notice.

#### Parameters

- **source** – absolute path to the source file on the DSO filesystem

- **destination** – absolute path to the destination file on the DSO filesystem

**Raises**

[RT01024Error](#) – if the operation did not complete

**get\_acquire\_length()** → float

Gets the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]

**Returns**

the time for one acquisition. Range: 250e-12 ... 500 [s]

**get\_channel\_offset(channel: int)** → float

Gets the voltage offset of the indicated channel.

**Parameters**

**channel** – is the channel number (1..4)

**Returns**

channel offset voltage in V (value between -1 and 1)

**get\_channel\_position(channel: int)** → float

Gets the vertical position of the indicated channel.

**Parameters**

**channel** – is the channel number (1..4)

**Returns**

channel position in div (value between -5 and 5)

**get\_channel\_range(channel: int)** → float

Queries the channel range in V.

**Parameters**

**channel** – is the input channel (1..4)

**Returns**

channel range in V

**get\_channel\_scale(channel: int)** → float

Queries the channel scale in V/div.

**Parameters**

**channel** – is the input channel (1..4)

**Returns**

channel scale in V/div

**get\_channel\_state(channel: int)** → bool

Queries if the channel is active or not.

**Parameters**

**channel** – is the input channel (1..4)

**Returns**

True if active, else False

**get\_reference\_point()** → int

Gets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]

**Returns**

the reference in %

**get\_repetitions()** → int

Get the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- \*RST = 1

**Returns**

the number of waveforms to acquire

**get\_timestamps()** → List[float]

Gets the timestamps of all recorded frames in the history and returns them as a list of floats.

**Returns**

list of timestamps in [s]

**Raises**

**RTO1024Error** – if the timestamps are invalid

**list\_directory(path: str)** → List[Tuple[str, str, int]]

List the contents of a given directory on the oscilloscope filesystem.

**Parameters**

**path** – is the path to a folder

**Returns**

a list of filenames in the given folder

**load\_configuration(filename: str)** → *None*

Load current settings from a configuration file. The filename has to be specified without base directory and ‘.dfl’ extension.

**Information from the manual** *ReCaLL* calls up the instrument settings from an intermediate memory identified by the specified number. The instrument settings can be stored to this memory using the command \*SAV with the associated number. It also activates the instrument settings which are stored in a file and loaded using *MMEMemory:LOAD:STATE* .

**Parameters**

**filename** – is the name of the settings file without path and extension

**local\_display(state: bool)** → *None*

Enable or disable local display of the scope.

**Parameters**

**state** – is the desired local display state

**prepare\_ultra\_segmentation()** → *None*

Make ready for a new acquisition in ultra segmentation mode. This function does one acquisition without ultra segmentation to clear the history and prepare for a new measurement.

**read\_measurement(meas\_n: int, name: str)** → float

**Parameters**

- **meas\_n** – measurement number 1..8
- **name** – measurement name, for example “MAX”

**Returns**

measured value

**run\_continuous\_acquisition()** → *None*

Start acquiring continuously.

**run\_single\_acquisition()** → *None*

Start a single or Nx acquisition.

**save\_configuration(filename: str)** → *None*

Save the current oscilloscope settings to a file. The filename has to be specified without path and ‘.dfl’ extension, the file will be saved to the configured settings directory.

**Information from the manual** SAVe stores the current instrument settings under the specified number in an intermediate memory. The settings can be recalled using the command \*RCL with the associated number. To transfer the stored instrument settings to a file, use *MMEMory:STORe:STATe* .

**Parameters**

**filename** – is the name of the settings file without path and extension

**save\_waveform\_history(filename: str, channel: int, waveform: int = 1)** → *None*

Save the history of one channel and one waveform to a .bin file. This function is used after an acquisition using sequence trigger mode (with or without ultra segmentation) was performed.

**Parameters**

- **filename** – is the name (without extension) of the file
- **channel** – is the channel number
- **waveform** – is the waveform number (typically 1)

**Raises**

**RT01024Error** – if storing waveform times out

**set\_acquire\_length(timerange: float)** → *None*

Defines the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]
- \*RST = 0.5 [s]

**Parameters**

**timerange** – is the time for one acquisition. Range: 250e-12 ... 500 [s]

**set\_channel\_offset**(channel: int, offset: float) → *None*

Sets the voltage offset of the indicated channel.

- Range: Dependent on the channel scale and coupling [V]
- Increment: Minimum 0.001 [V], may be higher depending on the channel scale and coupling
- \*RST = 0

**Parameters**

- **channel** – is the channel number (1..4)
- **offset** – Offset voltage. Positive values move the waveform down, negative values move it up.

**set\_channel\_position**(channel: int, position: float) → *None*

Sets the vertical position of the indicated channel as a graphical value.

- Range: -5.0 ... 5.0 [div]
- Increment: 0.02
- \*RST = 0

**Parameters**

- **channel** – is the channel number (1..4)
- **position** – is the position. Positive values move the waveform up, negative values move it down.

**set\_channel\_range**(channel: int, v\_range: float) → *None*

Sets the voltage range across the 10 vertical divisions of the diagram. Use the command alternatively instead of set\_channel\_scale.

- Range for range: Depends on attenuation factors and coupling. With 1:1 probe and external attenuations and 50 input coupling, the range is 10 mV to 10 V. For 1 M input coupling, it is 10 mV to 100 V. If the probe and/or external attenuation is changed, multiply the range values by the attenuation factors.
- Increment: 0.01
- \*RST = 0.5

**Parameters**

- **channel** – is the channel number (1..4)
- **v\_range** – is the vertical range [V]

**set\_channel\_scale**(channel: int, scale: float) → *None*

Sets the vertical scale for the indicated channel. The scale value is given in volts per division.

- Range for scale: depends on attenuation factor and coupling. With 1:1 probe and external attenuations and 50 input coupling, the vertical scale (input sensitivity) is 1 mV/div to 1 V/div. For 1 M input coupling, it is 1 mV/div to 10 V/div. If the probe and/or external attenuation is changed, multiply the values by the attenuation factors to get the actual scale range.
- Increment: 1e-3
- \*RST = 0.05

See also: `set_channel_range`

#### Parameters

- **channel** – is the channel number (1..4)
- **scale** – is the vertical scaling [V/div]

**set\_channel\_state**(*channel: int, state: bool*) → *None*

Switches the channel signal on or off.

#### Parameters

- **channel** – is the input channel (1..4)
- **state** – is True for on, False for off

**set\_reference\_point**(*percentage: int*) → *None*

Sets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]
- \*RST = 50 [%]

#### Parameters

**percentage** – is the reference in %

**set\_repetitions**(*number: int*) → *None*

Set the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- \*RST = 1

#### Parameters

**number** – is the number of waveforms to acquire

**set\_trigger\_level**(*channel: int, level: float, event\_type: int = 1*) → *None*

Sets the trigger level for the specified event and source.

- Range: -10 to 10 V
- Increment: 1e-3 V
- \*RST = 0 V

#### Parameters

- **channel** – indicates the trigger source.
  - 1..4 = channel 1 to 4, available for all event types 1..3
  - 5 = external trigger input on the rear panel for analog signals, available for A-event type = 1
  - 6..9 = not available
- **level** – is the voltage for the trigger level in [V].

- **event\_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

**set\_trigger\_mode(mode: Union[str, TriggerModes]) → None**

Sets the trigger mode which determines the behavior of the instrument if no trigger occurs.

#### Parameters

**mode** – is either auto, normal, or freerun.

#### Raises

**RT01024Error** – if an invalid triggermode is selected

**set\_trigger\_source(channel: int, event\_type: int = 1) → None**

Set the trigger (Event A) source channel.

#### Parameters

- **channel** – is the channel number (1..4)

- **event\_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

**start() → None**

Start the RT01024 oscilloscope and bring it into a defined state and remote mode.

**stop() → None**

Stop the RT01024 oscilloscope, reset events and close communication. Brings back the device to a state where local operation is possible.

**stop\_acquisition() → None**

Stop any acquisition.

```
class RT01024Config(waveforms_path: str, settings_path: str, backup_path: str, spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2, command_timeout_seconds: Union[int, float] = 60, wait_sec_short_pause: Union[int, float] = 0.1, wait_sec_enable_history: Union[int, float] = 1, wait_sec_post_acquisition_start: Union[int, float] = 2)
```

Bases: *VisaDeviceConfig*, *\_RT01024ConfigDefaultsBase*, *\_RT01024ConfigBase*

Configdataclass for the RT01024 device.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

#### Returns

a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### Returns

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**exception RT01024Error**

Bases: *DeviceError*

**class RT01024VisaCommunication(configuration)**

Bases: *VisaCommunication*

Specialization of VisaCommunication for the RTO1024 oscilloscope

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class RT01024VisaCommunicationConfig(host: Union[str, IPv4Address, IPv6Address], interface\_type:**

*Union[str, InterfaceType] = InterfaceType.TCPPIP\_INSTR, board: int = 0, port: int = 5025, timeout: int = 5000, chunk\_size: int = 204800, open\_timeout: int = 1000, write\_termination: str = '\n', read\_termination: str = '\n', visa\_backend: str = ''*

Bases: *VisaCommunicationConfig*

Configuration dataclass for VisaCommunication with specifications for the RTO1024 device class.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**interface\_type: Union[str, InterfaceType] = 2**

Interface type of the VISA connection, being one of *InterfaceType*.

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**


a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

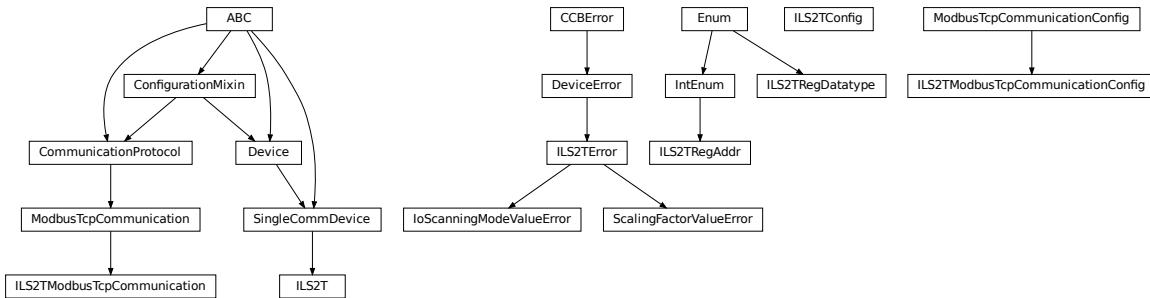**hvl\_ccb.dev.se\_ilst2t**

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

**class ILS2T(*com, dev\_config=None*)**

Bases: *SingleCommDevice*

Schneider Electric ILS2T stepper drive class.

**ACTION\_JOG\_VALUE = 0**

The single action value for *ILS2T.Mode.JOG*

**class ActionsPtp(*value*)**

Bases: *IntEnum*

Allowed actions in the point to point mode (*ILS2T.Mode.PTP*).

**ABSOLUTE\_POSITION = 0**

**RELATIVE\_POSITION\_MOTOR = 2**

**RELATIVE\_POSITION\_TARGET = 1**

**DEFAULT\_IO\_SCANNING\_CONTROL\_VALUES = {*'action': 2, 'continue\_after\_stop\_cu': 0, 'disable\_driver\_di': 0, 'enable\_driver\_en': 0, 'execute\_stop\_sh': 0, 'fault\_reset\_fr': 0, 'mode': 3, 'quick\_stop\_qs': 0, 'ref\_16': 1500, 'ref\_32': 0, 'reset\_stop\_ch': 0*}**

Default IO Scanning control mode values

**class Mode(*value*)**

Bases: *IntEnum*

ILS2T device modes

**JOG = 1**

**PTP = 3**

**class Ref16Jog(*value*)**

Bases: *Flag*

Allowed values for ILS2T ref\_16 register (the shown values are the integer representation of the bits), all in Jog mode = 1

**FAST** = 4

**NEG** = 2

**NEG\_FAST** = 6

**NONE** = 0

**POS** = 1

**POS\_FAST** = 5

#### **RegAddr**

Modbus Register Adresses

alias of [\*ILS2TRegAddr\*](#)

#### **RegDatatype**

Modbus Register Datatypes

alias of [\*ILS2TRegDatatype\*](#)

#### **class State(*value*)**

Bases: IntEnum

State machine status values

**ON** = 6

**QUICKSTOP** = 7

**READY** = 4

#### **static config\_cls()**

Return the default configdataclass class.

##### **Returns**

a reference to the default configdataclass class

#### **static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

##### **Returns**

the type of the standard communication protocol for this device

#### **disable(log\_warn: bool = True, wait\_sec\_max: Optional[int] = None) → bool**

Disable the driver of the stepper motor and enable the brake.

Note: the driver cannot be disabled if the motor is still running.

##### **Parameters**

- **log\_warn** – if log a warning in case the motor cannot be disabled.
- **wait\_sec\_max** – maximal wait time for the motor to stop running and to disable it; by default, with *None*, use a config value

##### **Returns**

*True* if disable request could and was sent, *False* otherwise.

**do\_iосcanning\_write**(\*\**kwargs*: int) → *None*

Perform a write operation using IO Scanning mode.

**Parameters**

**kwargs** – Keyword-argument list with options to send, remaining are taken from the defaults.

**enable()** → *None*

Enable the driver of the stepper motor and disable the brake.

**execute\_absolute\_position**(*position*: int) → bool

Execute a absolute position change, i.e. enable motor, perform absolute position change, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

**Parameters**

**position** – absolute position of motor in user defined steps.

**Returns**

*True* if actual position is as expected, *False* otherwise.

**execute\_relative\_step**(*steps*: int) → bool

Execute a relative step, i.e. enable motor, perform relative steps, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

**Parameters**

**steps** – Number of steps.

**Returns**

*True* if actual position is as expected, *False* otherwise.

**get\_dc\_volt()** → float

Read the DC supply voltage of the motor.

**Returns**

DC input voltage.

**get\_error\_code()** → Dict[int, Dict[str, Any]]

Read all messages in fault memory. Will read the full error message and return the decoded values. At the end the fault memory of the motor will be deleted. In addition, reset\_error is called to re-enable the motor for operation.

**Returns**

Dictionary with all information

**get\_position()** → int

Read the position of the drive and store into status.

**Returns**

Position step value

**get\_status()** → Dict[str, int]

Perform an IO Scanning read and return the status of the motor.

**Returns**

dict with status information.

**get\_temperature()** → int

Read the temperature of the motor.

**Returns**

Temperature in degrees Celsius.

**jog\_run(direction: bool = True, fast: bool = False)** → *None*

Slowly turn the motor in positive direction.

**jog\_stop()** → *None*

Stop turning the motor in Jog mode.

**quickstop()** → *None*

Stops the motor with high deceleration rate and falls into error state. Reset with *reset\_error* to recover into normal state.

**reset\_error()** → *None*

Resets the motor into normal state after quick stop or another error occurred.

**set\_jog\_speed(slow: int = 60, fast: int = 180)** → *None*

Set the speed for jog mode. Default values correspond to startup values of the motor.

**Parameters**

- **slow** – RPM for slow jog mode.
- **fast** – RPM for fast jog mode.

**set\_max\_acceleration(rpm\_minute: int)** → *None*

Set the maximum acceleration of the motor.

**Parameters**

**rpm\_minute** – revolution per minute per minute

**set\_max\_deceleration(rpm\_minute: int)** → *None*

Set the maximum deceleration of the motor.

**Parameters**

**rpm\_minute** – revolution per minute per minute

**set\_max\_rpm(rpm: int)** → *None*

Set the maximum RPM.

**Parameters**

**rpm** – revolution per minute ( 0 < rpm <= RPM\_MAX)

**Raises**

**ILS2TError** – if RPM is out of range

**set\_ramp\_type(ramp\_type: int = -1)** → *None*

**Set the ramp type. There are two options available:**

0: linear ramp -1: motor optimized ramp

**Parameters**

**ramp\_type** – 0: linear ramp | -1: motor optimized ramp

**start()** → *None*

Start this device.

**stop()** → *None*

Stop this device. Disables the motor (applies brake), disables access and closes the communication protocol.

**user\_steps(steps: int = 16384, revolutions: int = 1)** → *None*

Define steps per revolution. Default is 16384 steps per revolution. Maximum precision is 32768 steps per revolution.

**Parameters**

- **steps** – number of steps in *revolutions*.
- **revolutions** – number of revolutions corresponding to *steps*.

**write\_absolute\_position(position: int)** → *None*

Write instruction to turn the motor until it reaches the absolute position. This function does not enable or disable the motor automatically.

**Parameters**

**position** – absolute position of motor in user defined steps.

**write\_relative\_step(steps: int)** → *None*

Write instruction to turn the motor the relative amount of steps. This function does not enable or disable the motor automatically.

**Parameters**

**steps** – Number of steps to turn the motor.

```
class ILS2TConfig(rpm_max_init: Integral = 1500, wait_sec_post_enable: Union[int, float] = 1,
                    wait_sec_max_disable: Union[int, float] = 10, wait_sec_post_cannot_disable: Union[int,
                    float] = 1, wait_sec_post_relative_step: Union[int, float] = 2,
                    wait_sec_post_absolute_position: Union[int, float] = 2)
```

Bases: `object`

Configuration for the ILS2T stepper motor device.

**clean\_values()****force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True****classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**rpm\_max\_init: Integral = 1500**

initial maximum RPM for the motor, can be set up to 3000 RPM. The user is allowed to set a new max RPM at runtime using [ILS2T.set\\_max\\_rpm\(\)](#), but the value must never exceed this configuration setting.

**wait\_sec\_max\_disable: Union[int, float] = 10**

**wait\_sec\_post\_absolute\_position: Union[int, float] = 2**

**wait\_sec\_post\_CANNOT\_DISABLE: Union[int, float] = 1**

**wait\_sec\_post\_enable: Union[int, float] = 1**

**wait\_sec\_post\_relative\_step: Union[int, float] = 2**

**exception ILS2TError**

Bases: [DeviceError](#)

Error to indicate problems with the SE ILS2T stepper motor.

**class ILS2TModbusTcpCommunication(configuration)**

Bases: [ModbusTcpCommunication](#)

Specific implementation of Modbus/TCP for the Schneider Electric ILS2T stepper motor.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class ILS2TModbusTcpCommunicationConfig(host: Union[str, IPv4Address, IPv6Address], unit: int = 255, port: int = 502)**

Bases: [ModbusTcpCommunicationConfig](#)

Configuration dataclass for Modbus/TCP communication specific for the Schneider Electric ILS2T stepper motor.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**unit: int = 255**

The unit has to be 255 such that IO scanning mode works.

**class ILS2TRegAddr(value)**

Bases: IntEnum

Modbus Register Adresses for Schneider Electric ILS2T stepper drive.

**ACCESS\_ENABLE = 282**

**FLT\_INFO = 15362**

**FLT\_MEM\_DEL = 15112**

**FLT\_MEM\_RESET = 15114**

**IO\_SCANNING = 6922**

**JOGN\_FAST = 10506**

**JOGN\_SLOW = 10504**

**POSITION = 7706**

**RAMP\_ACC = 1556**

**RAMP\_DECEL = 1558**

**RAMP\_N\_MAX = 1554**

**RAMP\_TYPE = 1574**

**SCALE = 1550**

**TEMP = 7200**

**VOLT = 7198**

```
class ILS2TRegDatatype(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: `Enum`

Modbus Register Datatypes for Schneider Electric ILS2T stepper drive.

From the manual of the drive:

datatype	byte	min	max
INT8	1 Byte	-128	127
UINT8	1 Byte	0	255
INT16	2 Byte	-32_768	32_767
UINT16	2 Byte	0	65_535
INT32	4 Byte	-2_147_483_648	2_147_483_647
UINT32	4 Byte	0	4_294_967_295
BITS	just 32bits	N/A	N/A

`INT32 = (-2147483648, 2147483647)`

`is_in_range(value: int) → bool`

**exception IoScanningModeValueError**

Bases: `ILS2TError`


Error to indicate that the selected IO scanning mode is invalid.

**exception ScalingFactorValueError**

Bases: `ILS2TError`

Error to indicate that a scaling factor value is invalid.

## **hvl\_ccb.dev.sst\_luminox**

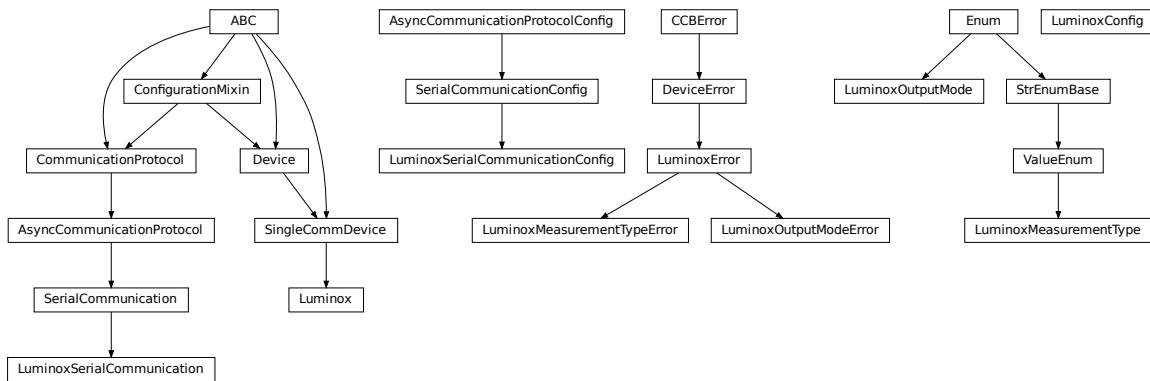

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.

Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

### `class Luminox(com, dev_config=None)`

Bases: `SingleCommDevice`

Luminox oxygen sensor device class.

#### `activate_output(mode: LuminoxOutputMode) → None`

activate the selected output mode of the Luminox Sensor. :param mode: polling or streaming

#### `static config_cls()`

Return the default configdataclass class.

##### **Returns**

a reference to the default configdataclass class

#### `static default_com_cls()`

Get the class for the default communication protocol used with this device.

##### **Returns**

the type of the standard communication protocol for this device

#### `query_polling(measurement: Union[str, LuminoxMeasurementType]) → Union[Dict[Union[str, LuminoxMeasurementType], Union[float, int, str]], float, int, str]`

Query a value or values of Luminox measurements in the polling mode, according to a given measurement type.

##### **Parameters**

`measurement` – type of measurement

##### **Returns**

value of requested measurement

##### **Raises**

- `ValueError` – when a wrong key for LuminoxMeasurementType is provided
- `LuminoxOutputModeError` – when polling mode is not activated
- `LuminoxMeasurementTypeError` – when expected measurement value is not read

#### `read_streaming() → Dict[Union[str, LuminoxMeasurementType], Union[float, int, str]]`

Read values of Luminox in the streaming mode. Convert the single string into separate values.

##### **Returns**

dictionary with `LuminoxMeasurementType.all_measurements_types()` keys and accordingly type-parsed values.

##### **Raises**

- `LuminoxOutputModeError` – when streaming mode is not activated
- `LuminoxMeasurementTypeError` – when any of expected measurement values is not read

#### `start() → None`

Start this device. Opens the communication protocol.

#### `stop() → None`

Stop the device. Closes also the communication protocol.

```
class LuminoxConfig(wait_sec_post_activate: Union[int, float] = 0.5, wait_sec_trials_activate: Union[int, float] = 0.1, nr_trials_activate: int = 5)
```

Bases: `object`

Configuration for the SST Luminox oxygen sensor.

`clean_values()`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

`is_configdataclass = True`

`classmethod keys() → Sequence[str]`

Returns a list of all configdataclass fields key-names.

#### Returns

a list of strings containing all keys.

`nr_trials_activate: int = 5`

`classmethod optional_defaults() → Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

#### Returns

a list of strings containing all optional keys.

`classmethod required_keys() → Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

#### Returns

a list of strings containing all required keys.

`wait_sec_post_activate: Union[int, float] = 0.5`

`wait_sec_trials_activate: Union[int, float] = 0.1`

`exception LuminoxError`

Bases: `DeviceError`

General Error for Luminox Device.

```
class LuminoxMeasurementType(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: `ValueEnum`

Measurement types for `LuminoxOutputMode.polling`.

The `all_measurements` type will read values for the actual measurement types as given in `LuminoxOutputMode.all_measurements_types()`; it parses multiple single values using regexp's for other measurement types, therefore, no regexp is defined for this measurement type.

```

all_measurements = 'A'

classmethod all_measurements_types() → Tuple[LuminoxMeasurementType, ...]

A tuple of LuminoxMeasurementType enum instances which are actual measurements, i.e. not date of manufacture or software revision.

barometric_pressure = 'P'

property command: str

date_of_manufacture = '# 0'

parse_read_measurement_value(read_txt: str) → Union[Dict[Union[str, LuminoxMeasurementType], Union[float, int, str]], float, int, str]

partial_pressure_o2 = '0'

percent_o2 = '%'

sensor_status = 'e'

serial_number = '# 1'

software_revision = '# 2'

temperature_sensor = 'T'

```

### **LuminoxMeasurementTypeDict**

A typing hint for a dictionary holding *LuminoxMeasurementType* values. Keys are allowed as strings because *LuminoxMeasurementType* is of a *StrEnumBase* type.

alias of Dict[Union[str, LuminoxMeasurementType], Union[float, int, str]]

### **exception LuminoxMeasurementTypeError**

Bases: *LuminoxError*

Wrong measurement type for requested data

### **LuminoxMeasurementTypeValue**

A typing hint for all possible *LuminoxMeasurementType* values as read in either streaming mode or in a polling mode with *LuminoxMeasurementType.all\_measurements*.

Beware: has to be manually kept in sync with *LuminoxMeasurementType* instances *cast\_type* attribute values.

alias of Union[float, int, str]

### **class LuminoxOutputMode(value)**

Bases: *Enum*

output mode.

**polling** = 1

**streaming** = 0

### **exception LuminoxOutputModeError**

Bases: *LuminoxError*

Wrong output mode for requested data

**class LuminoxSerialCommunication**(*configuration*)

Bases: *SerialCommunication*

Specific communication protocol implementation for the SST Luminox oxygen sensor. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**class LuminoxSerialCommunicationConfig**(*terminator: bytes = b'\r\n'*, *encoding: str = 'utf-8'*,

*encoding\_error\_handling: str = 'strict'*,

*wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5*,

*default\_n\_attempts\_read\_text\_nonempty: int = 10*, *port: Union[str, NoneType] = None*,

*baudrate: int = 9600*, *parity: Union[SerialCommunicationParity, NoneType] = <SerialCommunicationParity.NONE: 'N'>*,

*stopbits: Union[int, SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>*,

*bytesize: Union[int, SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>*,

*timeout: Union[int, float] = 3*)

Bases: *SerialCommunicationConfig*

**baudrate: int = 9600**

Baudrate for SST Luminox is 9600 baud

**bytesize: Union[int, SerialCommunicationBytesize] = 8**

One byte is eight bits long

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**parity: Union[str, SerialCommunicationParity] = 'N'**

SST Luminox does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**stopbits: Union[int, SerialCommunicationStopbits] = 1**

SST Luminox does use one stop bit

**terminator: bytes = b'\r\n'**

The terminator is CR LF

**timeout: Union[int, float] = 3**

use 3 seconds timeout as default

## hvl\_ccb.dev.utils

### Poller

**class Poller(spoll\_handler: Callable, polling\_delay\_sec: Union[int, float] = 0, polling\_interval\_sec: Union[int, float] = 1, polling\_timeout\_sec: Optional[Union[int, float]] = None)**

Bases: object

Poller class wrapping *concurrent.futures.ThreadPoolExecutor* which enables passing of results and errors out of the polling thread.

**is\_polling()** → bool

Check if device status is being polled.

**Returns**

*True* when polling thread is set and alive

**start\_polling()** → bool

Start polling.

**Returns**

*True* if was not polling before, *False* otherwise

**stop\_polling()** → bool

Stop polling.

Wait for until polling function returns a result as well as any exception / error that might have been raised within a thread.

**Returns**

*True* if was polling before, *False* otherwise, and last result of the polling function call.

**Raises**

polling function exceptions

**wait\_for\_polling\_result()**


Wait for until polling function returns a result as well as any exception / error that might have been raised within a thread.

**Returns**

polling function result

**Raises**

polling function errors

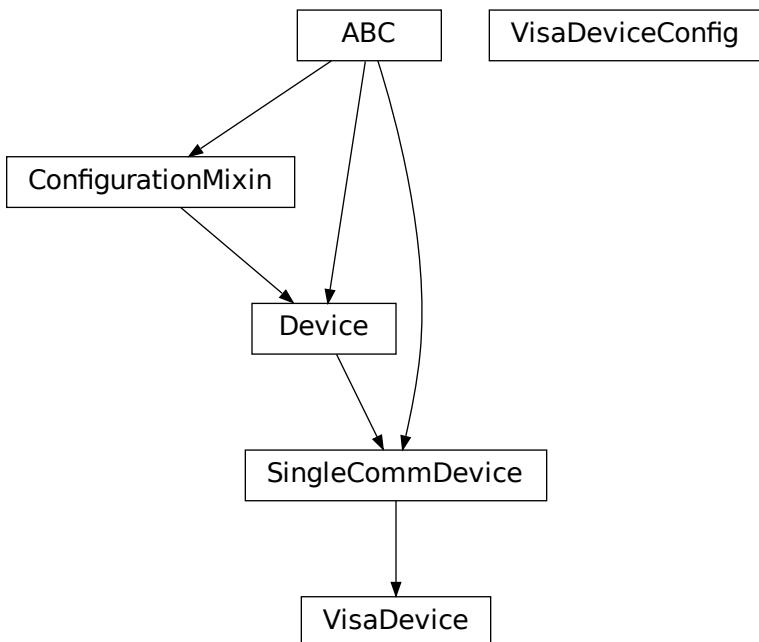**hvl\_ccb.dev.visa**

```
class VisaDevice(com: Union[VisaCommunication, VisaCommunicationConfig, dict], dev_config: Optional[Union[VisaDeviceConfig, dict]] = None)
```

Bases: `SingleCommDevice`

Device communicating over the VISA protocol using `VisaCommunication`.

**static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**static default\_com\_cls() → Type[VisaCommunication]**

Return the default communication protocol for this device type, which is `VisaCommunication`.

**Returns**

the VisaCommunication class

**get\_error\_queue() → str**

Read out error queue and logs the error.

**Returns**

Error string

**get\_identification() → str**

Queries “\*IDN?” and returns the identification string of the connected device.

**Returns**

the identification string of the connected device

**reset() → None**

Send “\*RST” and “\*CLS” to the device. Typically sets a defined state.

**spoll\_handler()**

Reads the status byte and decodes it. The status byte STB is defined in IEEE 488.2. It provides a rough overview of the instrument status.

**Returns****start() → None**

Start the VisaDevice. Sets up the status poller and starts it.

**Returns****stop() → None**

Stop the VisaDevice. Stops the polling thread and closes the communication protocol.

**Returns****wait\_operation\_complete(timeout: Optional[float] = None) → bool**

Waits for a operation complete event. Returns after timeout [s] has expired or the operation complete event has been caught.

**Parameters**

**timeout** – Time in seconds to wait for the event; *None* for no timeout.

**Returns**

True, if OPC event is caught, False if timeout expired

**class VisaDeviceConfig(spoll\_interval: Union[int, float] = 0.5, spoll\_start\_delay: Union[int, float] = 2)**

Bases: `_VisaDeviceConfigDefaultsBase`, `_VisaDeviceConfigBase`

Configdataclass for a VISA device.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

## Module contents

Devices subpackage.

### hvl\_ccb.utils

#### Subpackages

##### hvl\_ccb.utils.conversion

#### Submodules

##### hvl\_ccb.utils.conversion.map\_range

MapBitAsymRange

MapBitSymRange

MapRanges

**class MapBitAsymRange**(value: ~typing.Union[int, float], bit: int, dtype\_1: ~typing.Union[~numpy.dtype, None, type, ~numpy.typing.\_dtype\_like.\_SupportsDType[~numpy.dtype]], str, ~typing.Tuple[~typing.Any, int], ~typing.Tuple[~typing.Any, ~typing.Union[~typing\_extensions.SupportsIndex, ~typing.Sequence[~typing\_extensions.SupportsIndex]]], ~typing.List[~typing.Any], ~numpy.typing.\_dtype\_like.\_DTypeDict, ~typing.Tuple[~typing.Any, ~typing.Any]] = <class 'float'>, logger=None)

Bases: `_MapBitRange`

Class to convert an asymmetric arbitrary range (0 to value) to a bit-range (0 to  $2^{**\text{bit}} - 1$ ).

```
class MapBitSymRange(value: ~typing.Union[int, float], bit: int, dtype_1: ~typing.Union[~numpy.dtype, None,
    type, ~numpy.typing._dtype_like._SupportsDType[~numpy.dtype], str,
    ~typing.Tuple[~typing.Any, int], ~typing.Tuple[~typing.Any,
    ~typing.Union[~typing_extensions.SupportsIndex,
    ~typing.Sequence[~typing_extensions.SupportsIndex]]], ~typing.List[~typing.Any],
    ~numpy.typing._dtype_like._DTypeDict, ~typing.Tuple[~typing.Any, ~typing.Any]] = <class 'float'>, logger=None)
```

Bases: `_MapBitRange`

Class to convert a symmetric arbitrary range (-value to value) to a bit-range (0 to  $2^{**\text{bit}} - 1$ ).


```
class MapRanges(range_1: Tuple[Union[int, float], Union[int, float]], range_2: Tuple[Union[int, float],
    Union[int, float]], dtype_1: Union[dtype, None, type, _SupportsDType[dtype], str, Tuple[Any,
    int], Tuple[Any, Union[SupportsIndex, Sequence[SupportsIndex]]], List[Any], _DTypeDict,
    Tuple[Any, Any]], dtype_2: Union[dtype, None, type, _SupportsDType[dtype], str, Tuple[Any,
    int], Tuple[Any, Union[SupportsIndex, Sequence[SupportsIndex]]], List[Any], _DTypeDict,
    Tuple[Any, Any]], logger=None)
```

Bases: `object`

`convert_to_range1(value, **kwargs)`

`convert_to_range2(value, **kwargs)`

## `hvl_ccb.utils.conversion.sensor`

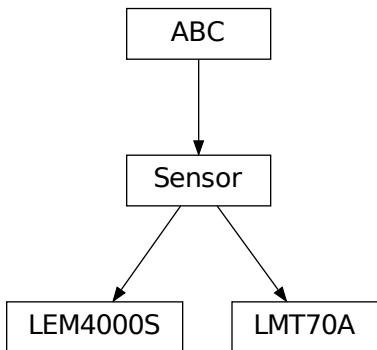

Sensors that are used by the devices implemented in the CCB

```
class LEM4000S
```

Bases: `Sensor`

`CONVERSION`

`calibration_factor`

`convert(value, **kwargs)`

`shunt`

```
class LMT70A
```

Bases: *Sensor*

Converts the output voltage (V) to the measured temperature (default °C) when using a TI Precision Analog Temperature Sensor LMT70(A)

LUT

```
convert(value, **kwargs)
```


```
temperature_unit
```

```
class Sensor
```

Bases: ABC

```
abstract convert(value, **kwargs)
```

[hvl\\_ccb.utils.conversion.unit](#)

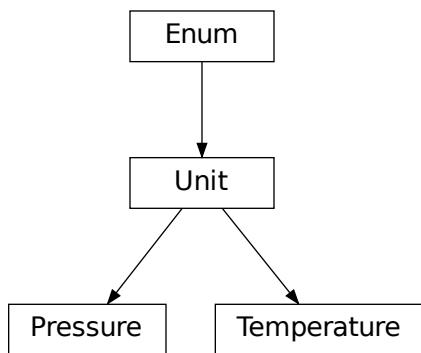

Unit conversion, within in the same group of units, for example Kelvin <-> Celsius

```
class Pressure(value)
```

Bases: *Unit*

An enumeration.

```
ATM = 'atm'
```

```
ATMOSPHERE = 'atm'
```

```
BAR = 'bar'
```

```
MILLIMETER_MERCURY = 'mmHg'
```

```
MMHG = 'mmHg'
```

```
PA = 'Pa'
```

```
PASCAL = 'Pa'
```

```

POUNDS_PER_SQUARE_INCH = 'psi'
PSI = 'psi'
TORR = 'torr'

class Temperature(value)
    Bases: Unit
    An enumeration.


    C = 'C'
    CELSIUS = 'C'
    F = 'F'
    FAHRENHEIT = 'F'
    K = 'K'
    KELVIN = 'K'

class Unit(value)
    Bases: Enum
    An enumeration.

    abstract classmethod convert(value, **kwargs)

```

## [hvl\\_ccb.utils.conversion.utils](#)

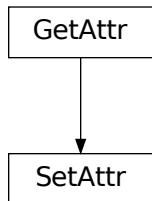

```

class GetAttr(default, name)
    Bases: object

class SetAttr(default, name, limits, absolut=False, dtype=(<class 'int'>, <class 'float'>), validator=None)
    Bases: GetAttr


convert_value_to_str(*value: ndarray[Any, dtype[ScalarType]]) → List[Union[str, List[str]]]

```

Converts two sets of values to strings. This is necessary because a 0-dim array needs different treatment than a 1-dim array :param value: array of values either 0-dim or 1-dim :return:converted se

**preserve\_type(func)**

This wrapper preserves the first order type of the input. Upto now the type of the data stored in a list, tuple, array or dict is not preserved. Integer will be converted to float!

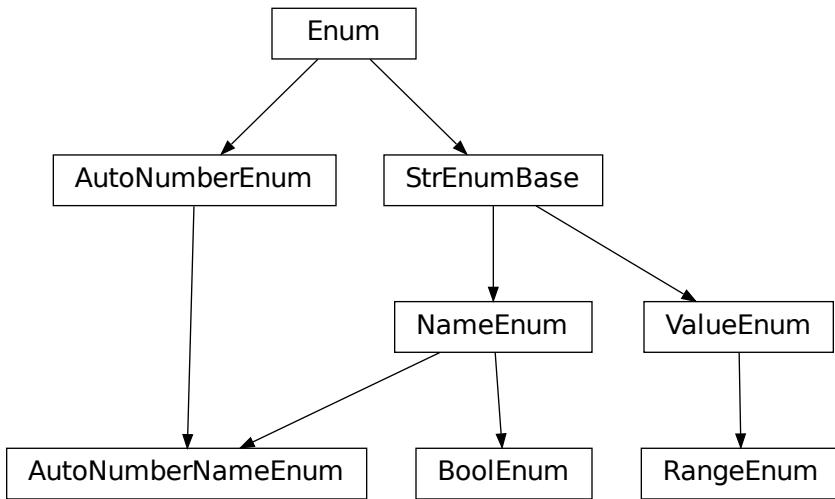**Module contents****Submodules****hvl\_ccb.utils.enum**

```
class AutoNumberNameEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: [NameEnum](#), [AutoNumberEnum](#)

Auto-numbered enum with names used as string representation, and with lookup and equality based on this representation.

```
class BoolEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: [NameEnum](#)

**BoolEnum inherits from NameEnum and the type of the first value is**  
enforced to be ‘boolean’. For bool()-operation the `__bool__` is redefined here.

```
class NameEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1, boundary=None)
```

Bases: [StrEnumBase](#)

Enum with names used as string representation, and with lookup and equality based on this representation. The lookup is implemented in `StrEnumBase` with the `_missing_value_` method. The equality is also defined at this place (`__eq__`).

Use-case:

```
class E(NameEnum):
    a = 2
    b = 4

E.a == "a"
E.a != 2
E.a != "2"
```

The access would be normally with E["a"], but E("a") works also. Therefore, E["a"] == E("a")

Attention: to avoid errors, best use together with *unique* enum decorator.

```
class RangeEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                boundary=None)
```

Bases: float, [ValueEnum](#)

Range enumeration inherit from ValueEnum, find suitable voltage/current/resistance input range for devices such as multimeter and oscilloscope

```
abstract classmethod unit() → str
```

Returns the Unit of the values in the enumeration. :return: the unit of the values in the enumeration in string format

```
class StrEnumBase(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                  boundary=None)
```

Bases: Enum

String representation-based equality and lookup.

```
class ValueEnum(value=<no_arg>, names=None, module=None, qualname=None, type=None, start=1,
                 boundary=None)
```

Bases: [StrEnumBase](#)

Enum with string representation of values used as string representation, and with lookup and equality based on this representation. Values do not need to be of type 'str', but they need to have a str-representation to enable this feature. The lookup is implemented in StrEnumBase with the `_missing_value_` method. The equality is also defined at this place (`__eq__`).

Use-case:

```
class E(ValueEnum):
    ONE = 1

E.ONE == "1"
E.ONE != 1
E.ONE != "ONE"
```

The access would be normally with E(1), but E("1") works also. Therefore, E(1) == E("1")

Attention: to avoid errors, best use together with *unique* enum decorator.

## **hvl\_ccb.utils.typing**

Additional Python typing module utilities

### **ConvertibleTypes**

Typing hint for data type that can be used in conversion

alias of Union[int, float, List[Union[int, float]], Tuple[Union[int, float], ...], Dict[str, Union[int, float]], ndarray]

### **Number**

Typing hint auxiliary for a Python base number types: *int* or *float*.

alias of Union[int, float]

### **check\_generic\_type(value, type\_, name='instance')**

Check if *value* is of a generic type *type\_*. Raises *TypeError* if it's not.

#### **Parameters**

- **name** – name to report in case of an error
- **value** – value to check
- **type** – generic type to check against

### **is\_generic\_type\_hint(type\_)**

Check if class is a generic type, for example *Union[int, float]*, *List[int]*, or *List*.

#### **Parameters**

**type** – type to check

## **hvl\_ccb.utils.validation**

### **validate\_and\_resolve\_host(host: Union[str, IPv4Address, IPv6Address], logger: Optional[Logger] = None) → str**

### **validate\_bool(x\_name: str, x: object, logger: Optional[Logger] = None) → None**

Validate if given input *x* is a *bool*.

#### **Parameters**

- **x\_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as boolean
- **logger** – logger of the calling submodule

#### **Raises**

**TypeError** – when the validated input does not have boolean type

### **validate\_number(x\_name: str, x: object, limits: ~typing.Optional[~typing.Tuple] = (None, None), number\_type: ~typing.Union[~typing.Type[~typing.Union[int, float]], ~typing.Tuple[~typing.Type[~typing.Union[int, float]], ...]] = (<class 'int'>, <class 'float'>), logger: ~typing.Optional[~logging.Logger] = None) → None**

Validate if given input *x* is a number of given *number\_type* type, with value between given *limits[0]* and *limits[1]* (inclusive), if not *None*. For array-like objects (npt.NDArray, List, Tuple, Dict) it is checked if all elements are within the limits and have the correct type.


#### **Parameters**

- **x\_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as number of given type within given range
- **logger** – logger of the calling submodule
- **limits** – [lower, upper] limit, with *None* denoting no limit: [-inf, +inf]
- **number\_type** – expected type or tuple of types of a number, by default (*int, float*)

**Raises**

- **TypeError** – when the validated input does not have expected type
- **ValueError** – when the validated input has correct number type but is not within given range

**validate\_tcp\_port**(*port: int, logger: Optional[Logger] = None*)

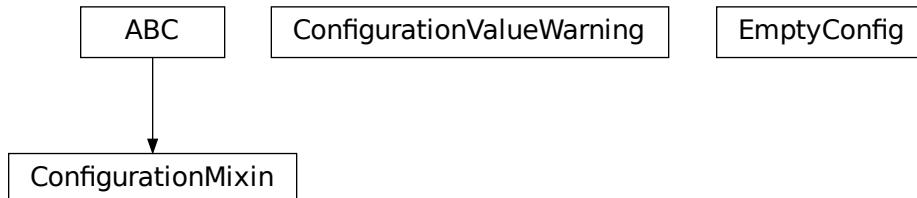**Module contents****4.1.2 Submodules****hvl\_ccb.configuration****Facilities**

providing classes for handling configuration for communication protocols and devices.

**class ConfigurationMixin(configuration)**

Bases: ABC

Mixin providing configuration to a class.

**property config**

ConfigDataclass property.

**Returns**

the configuration

**abstract static config\_cls()**

Return the default configdataclass class.

**Returns**

a reference to the default configdataclass class

**configuration\_save\_json**(*path*: str) → *None*

Save current configuration as JSON file.

**Parameters**

**path** – path to the JSON file.

**classmethod from\_json**(*filename*: str)

Instantiate communication protocol using configuration from a JSON file.

**Parameters**

**filename** – Path and filename to the JSON configuration

**exception ConfigurationValueWarning**

Bases: UserWarning

User warnings category for values of @configdataclass fields.

**class EmptyConfig**

Bases: object

Empty configuration dataclass.

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns**

a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns**

a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns**

a list of strings containing all required keys.

**configdataclass**(*direct\_decoration=None, frozen=True*)

Decorator to make a class a configdataclass. Types in these dataclasses are enforced. Implement a function `clean_values(self)` to do additional checking on value ranges etc.

It is possible to inherit from a configdataclass and re-decorate it with `@configdataclass`. In a subclass, default values can be added to existing fields. Note: adding additional non-default fields is prone to errors, since the order has to be respected through the whole chain (first non-default fields, only then default-fields).

**Parameters**


**frozen** – defaults to True. False allows to later change configuration values. Attention: if configdataclass is not frozen and a value is changed, typing is not enforced anymore!

**hvl\_ccb.error**

Introduce a common code base error for the CCB

**exception CCBError**

Bases: `Exception`

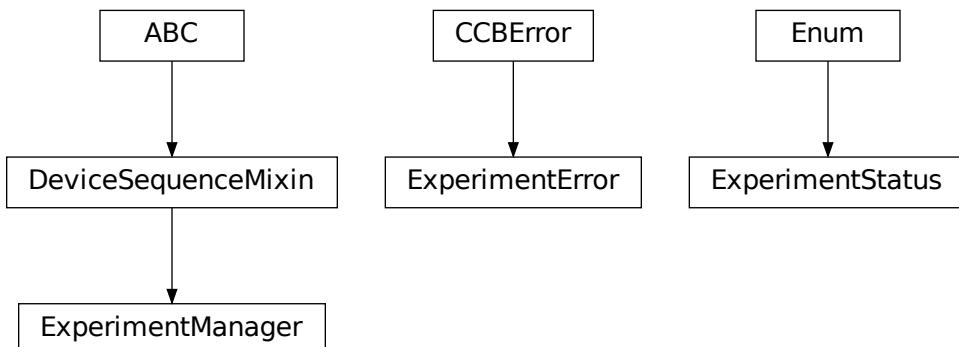**hvl\_ccb.experiment\_manager**

Main module containing the top level `ExperimentManager` class. Inherit from this class to implement your own experiment functionality in another project and it will help you start, stop and manage your devices.

**exception ExperimentError**

Bases: `CCBError`

Error to indicate that the current status of the experiment manager is on ERROR and thus no operations can be made until reset.

```
class ExperimentManager(*args, **kwargs)
```

Bases: *DeviceSequenceMixin*

Experiment Manager can start and stop communication protocols and devices. It provides methods to queue commands to devices and collect results.

**add\_device**(*name*: str, *device*: Device) → *None*

Add a new device to the manager. If the experiment is running, automatically start the device. If a device with this name already exists, raise an error.

**Parameters**

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

**Raises**

*ExperimentError* –

**devices\_failed\_start**: Dict[str, Device]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices\_failed\_stop*.

**devices\_failed\_stop**: Dict[str, Device]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices\_failed\_start*.

**finish**() → *None*

Stop experimental setup, stop all devices.

**is\_error**() → bool

Returns true, if the status of the experiment manager is *error*.

**Returns**

True if on error, false otherwise

**is\_finished**() → bool

Returns true, if the status of the experiment manager is *finished*.

**Returns**

True if finished, false otherwise

**is\_running**() → bool

Returns true, if the status of the experiment manager is *running*.

**Returns**

True if running, false otherwise

**run**() → *None*

Start experimental setup, start all devices.

**start**() → *None*

Alias for ExperimentManager.run()

**property status**: ExperimentStatus

Get experiment status.

**Returns**

experiment status enum code.

**stop()** → *None*

Alias for ExperimentManager.finish()

**class ExperimentStatus(*value*)**

Bases: `Enum`

Enumeration for the experiment status

**ERROR = 5**

**FINISHED = 4**

**FINISHING = 3**

**INITIALIZED = 0**

**INITIALIZING = -1**

**RUNNING = 2**

**STARTING = 1**

#### 4.1.3 Module contents

Top-level package for HVL Common Code Base.



## **CONTRIBUTING**

Contributions are welcome, and they are greatly appreciated! Every little bit helps, and credit will always be given.

You can contribute in many ways:

### **5.1 Types of Contributions**

#### **5.1.1 Report Bugs**

Report bugs at [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/issues](https://gitlab.com/ethz_hvl/hvl_ccb/issues).

If you are reporting a bug, please include:

- Your operating system name and version.
- Any details about your local setup that might be helpful in troubleshooting.
- Detailed steps to reproduce the bug.

#### **5.1.2 Fix Bugs**

Look through the GitLab issues for bugs. Anything tagged with “bug” and “help wanted” is open to whoever wants to implement it.

#### **5.1.3 Implement Features**

Look through the GitLab issues for features. Anything tagged with “enhancement” and “help wanted” is open to whoever wants to implement it.

#### **5.1.4 Write Documentation**

HVL Common Code Base could always use more documentation, whether as part of the official HVL Common Code Base docs, in docstrings, or even on the web in blog posts, articles, and such.

### 5.1.5 Submit Feedback

The best way to send feedback is to file an issue at [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/issues](https://gitlab.com/ethz_hvl/hvl_ccb/issues).

If you are proposing a feature:

- Explain in detail how it would work.
- Keep the scope as narrow as possible, to make it easier to implement.
- Remember that this is a volunteer-driven project, and that contributions are welcome :)

## 5.2 Get Started!

Ready to contribute? Here's how to set up *hvl\_ccb* for local development.

1. Clone *hvl\_ccb* repo from GitLab.

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

2. Install your local copy into a virtualenv. Assuming you have `virtualenvwrapper` installed, this is how you set up your fork for local development:

```
$ mkvirtualenv hvl_ccb
$ cd hvl_ccb/
$ pip install -e .[all]
$ pip install -r requirements_dev.txt
```

3. Create a branch for local development:

```
$ git checkout -b name-of-your-bugfix-or-feature
```

Now you can make your changes locally.

4. When you're done making changes, check that your changes pass flake8, mypy and the tests, including testing other Python versions with tox:

```
$ flake8 hvl_ccb tests
$ mypy --show-error-codes hvl_ccb
$ python setup.py test or py.test
$ tox
```

You can also use the provided make-like shell script to run flake8 and tests:

```
$ ./make.sh style
$ ./make.sh type
$ ./make.sh test
```

5. As we want to maintain a high quality of coding style we use *black*. This style is checked with the pipelines on gitlab.com. Ensure that your commits include only properly formatted code. One way to comply is to install and use *pre-commit*. This package includes the necessary configuration.

6. Commit your changes and push your branch to GitLab:

```
$ git add .
$ git commit -m "Your detailed description of your changes."
$ git push origin name-of-your-bugfix-or-feature
```

7. Submit a merge request through the GitLab website.

## 5.3 Merge Request Guidelines

Before you submit a merge request, check that it meets these guidelines:

1. The merge request should include tests.
2. If the merge request adds functionality, the docs should be updated. Put your new functionality into a function with a docstring, and add the feature to the list in README.rst.
3. The merge request should work for Python 3.7 to 3.10. Check [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/merge\\_requests](https://gitlab.com/ethz_hvl/hvl_ccb/merge_requests) and make sure that the tests pass for all supported Python versions.

## 5.4 Tips

- To run tests from a single file:

```
$ py.test tests/test_hvl_ccb.py
```

or a single test function:

```
$ py.test tests/test_hvl_ccb.py::test_command_line_interface
```

- If your tests are slow, profile them using the pytest-profiling plugin:

```
$ py.test tests/test_hvl_ccb.py --profile
```

or for a graphical overview (you need a SVG image viewer):

```
$ py.test tests/test_hvl_ccb.py --profile-svg
$ open prof/combined.svg
```

- To add dependency, edit appropriate \*requirements variable in the setup.py file and re-run:

```
$ python setup.py develop
```

- To generate a PDF version of the Sphinx documentation instead of HTML use:

```
$ rm -rf docs/hvl_ccb.rst docs/modules.rst docs/_build && sphinx-apidoc -o docs/hvl_ccb && python -msphinx -M latexpdf docs/ docs/_build
```

This command can also be run through the make-like shell script:

```
$ ./make.sh docs-pdf
```

This requires a local installation of a LaTeX distribution, e.g. MikTeX.

## 5.5 Deploying

A reminder for the maintainers on how to deploy.

Make sure all your changes are committed and that all relevant MR are merged. Then switch to `devel`, update it and create `release-N.M.K` branch:

```
$ git switch devel  
$ git pull  
$ git checkout -b release-N.M.K
```

- Update or create entry in `HISTORY.rst` (commit message: Update `HISTORY.rst`: release N.M.K).
- Update, if applicable, `AUTHORS.rst` (commit message: Update `AUTHORS.rst`: release N.M.K)
- Update features tables in `README.rst` file (commit message: Update `README.rst`: release N.M.K)
- Update API docs (commit message: Update API-docs: release N.M.K)

```
$ ./make.sh docs # windows  
$ make docs # unix-based-os
```

Commit all of the above, except for

- `docs/hvl_ccb.dev.picotech_pt104.rst`
- `docs/hvl_ccb.dev.tiepie.base.rst`
- `docs/hvl_ccb.dev.tiepie.channel.rst`
- `docs/hvl_ccb.dev.tiepie.device.rst`
- `docs/hvl_ccb.dev.tiepie.generator.rst`
- `docs/hvl_ccb.dev.tiepie.i2c.rst`
- `docs/hvl_ccb.dev.tiepie.oscilloscope.rst`
- `docs/hvl_ccb.dev.tiepie.utils.rst`.

Before you continue revert the changes in this file.

Then run:

```
$ bumpversion patch # possible: major / minor / patch  
$ git push --set-upstream origin release-N.M.K  
$ git push --tags
```

Go to <https://readthedocs.org/projects/hvlccb/builds/> and check if RTD docs build for the pushed tag passed.

Wait for the CI pipeline to finish successfully.

The two following commands are best executed in a WSL or Unix based OS. Run a release check:

```
$ make release-check
```

Finally, prepare and push a release:

```
$ make release
```

Merge the release branch into master and devel branches with `--no-ff` flag and delete the release branch:

```
$ git switch master
$ git merge --no-ff release-N.M.K
$ git push
$ git switch devel
$ git merge --no-ff release-N.M.K
$ git push
$ git push --delete origin release-N.M.K
$ git branch --delete release-N.M.K
```

After this you can/should clean your folder (with WSL/Unix command):

```
$ make clean
```

Finally, prepare GitLab release and cleanup the corresponding milestone:

1. go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/-/tags/](https://gitlab.com/ethz_hvl/hvl_ccb/-/tags/), select the latest release tag, press “Edit release notes” and add the release notes (copy a corresponding entry from `HISTORY.rst` file with formatting adjusted from ReStructuredText to Markdown); press “Save changes”;
2. go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/-/releases](https://gitlab.com/ethz_hvl/hvl_ccb/-/releases), select the latest release, press “Edit this release” and under “Milestones” select the corresponding milestone; press “Save changes”;
3. go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/-/milestones](https://gitlab.com/ethz_hvl/hvl_ccb/-/milestones), make sure that it is 100% complete (otherwise, create a next patch-level milestone and assign it to the ongoing Issues and Merge Requests therein); press “Close Milestone”.



**CREDITS**

## 6.1 Maintainers

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>
- Chi-Ching Hsu <hsu@eeh.ee.ethz.ch>

## 6.2 Authors

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- David Gruber <dev@davidgruber.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Alise Chachereau <chachereau@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- David Taylor <dtaylor@ethz.ch>
- Joseph Engelbrecht <engelbrecht@eeh.ee.ethz.ch>
- Chi-Ching Hsu <hsu@eeh.ee.ethz.ch>

## 6.3 Contributors

- Luca Nembrini <lucane@student.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>
- Raphael Faerber <raphael.faerber@eeh.ee.ethz.ch>
- Ruben Stadler <rstadler@student.ethz.ch>
- Hanut Vemulapalli <vemulapalli@eeh.ee.ethz.ch>



## HISTORY

### 7.1 0.12.1 (2022-10-31)

- Fix **numpy** version requirement problem
  - for Python 3.7: 1.21.6
  - for Python 3.8 and onwards: 1.23.4

### 7.2 0.12.0 (2022-10-17)

- Last release for Python 3.7 and 3.8
- **Repository maintenance**
  - update Labjack LJM software installer link in the pipeline
  - fix dependencies to the fixed version
  - fix `asyncua` to 0.9.95 and `pymodbus` to 2.5.3 (newer versions break the code)
  - fix PICube checker for slope as it is always positive

### 7.3 0.11.1 (2022-09-15)

- **Repository maintenance**
  - fix issue with `mypy` and Python 3.10.7
  - update code style to `black` 22.8.0
  - project configurations merged into `setup.cfg`
  - fix coverage indicator

## 7.4 0.11.0 (2022-06-22)

- New device: Fluke 884X Bench 6.5 Digit Precision Multimeter
- RangeEnum is a new enum for e.g. measurement ranges which also finds a suitable range object
- **smaller changes of device tiepie:**
  - introduce status method `is_measurement_running()` to check if the device is armed
  - introduce `stop_measurement()` to disarm the trigger of the device
  - fix bug with docs due to change of libtiepie
- NameEnum and inherited enums can only have unique entries

## 7.5 0.10.3 (2022-03-21)

- fix bug in the Labjack pulse feature that occurred when the start time was set to 0s
- new conversion utility to map two ranges on each other
- update CONTRIBUTING.RST
- update makefile and make.sh
- improve the mockup telnet test server

## 7.6 0.10.2 (2022-02-28)

- introduction of `black` as code formatter
- increase the required version of the package `aenum`
- remove device `supercube2015` - as it is no longer used
- remove unused package `openpyxl` requirement
- fix bug in highland logging
- improve handling for communication error with picotech

## 7.7 0.10.1 (2022-01-24)

- **several improvements and fixes for device cube:**
  - privatize `Alarms` and `AlarmsOverview`
  - fix list of cube alarms
  - improve docs
  - fix bugs with earthing sticks
  - fix bug in config dataclass of cube
- introduction of `BoolEnum`
- introduction of `RangeEnum`

- bumpversion -> bump2version

## 7.8 0.10.0 (2022-01-17)

- Reimplementation of the Cube (before known as Supercube)
- **new names:**
  - Supercube Typ B -> BaseCube
  - Supercube Typ A -> PICube (power inverter Cube)
- **new import:**
  - `from hvl_ccb.dev.supercube import SupercubeB` -> `from hvl_ccb.dev.cube import BaseCube`
- **new programming style:**
  - getter / setter methods -> properties
  - e.g. get: `cube.get_support_output(port=1, contact=1)` -> `cube.support_1.output_1`
  - e.g. set: `cube.get_support_output(port=1, contact=1, state=True)` -> `cube.support_1.output_1 = True`
- unify Exceptions of Cube
- implement Fast Switch-Off of Cube
- remove method `support_output_impulse`
- all active alarms can now be queried `cube.active_alarms()`
- alarms will now result in different logging levels depending on the seriousness of the alarm.
- introduction of limits for slope and safety limit for RedReady
- during the startup the CCB will update the time of the cube.
- verification of inputs
- polarity of DC voltage
- Switch from `python-opcua` to `opcua-asyncio` (former package is no longer maintained)

## 7.9 0.9.0 (2022-01-07)

- New device: Highland T560 digital delay and pulse generator over Telnet.
- **Rework of the Technix Capacitor Charger.**
  - Moved into a separate sub-package
  - NEW import over `import hvl_ccb.dev.technix as XXX`
  - Slightly adapted behaviour
- Add `validate_tcp_port` to validate port number.
- **Add `validate_and_resolve_host` to validate and resolve host names and IPs.**
  - Remove requirement IPy

- Add a unified CCB Exception schema for all devices and communication protocols.
- Add data conversion functions to README.
- Update CI and devel images from Debian 10 buster to Debian 11 bullseye.
- Fix typing due to numpy update.
- Fix incorrect overloading of `clean_values()` in classes of type `XCommunicationConfig`.

## 7.10 0.8.5 (2021-11-05)

- Added arbitrary waveform for TiePie signal generation, configurable via `dev.tiepie.generator.TiePieGeneratorConfig.waveform` property.
- In `utils.conversion_sensor`: improvements for class constants; removed SciPy dependency.
- Added Python 3.10 support.

## 7.11 0.8.4 (2021-10-22)

- `utils.validation.validate_number` extension to handle NumPy arrays and array-like objects.
- `utils.conversion_unit` utility classes handle correctly `NamedTuple` instances.
- `utils.conversion_sensor` and `utils.conversion_unit` code simplification (no `transfer_function_order` attribute) and cleanups.
- Fixed incorrect error logging in `configuration.configdataclass`.
- `comm.telnet.TelnetCommunication` tests fixes for local run errors.

## 7.12 0.8.3 (2021-09-27)

- New data conversion functions in `utils.conversion_sensor` and `utils.conversion_unit` modules. Note: to use these functions you must install `hvl_ccb` with extra requirement, either `hvl_ccb[conversion]` or `hvl_ccb[all]`.
- Improved documentation with respect to installation of external libraries.

## 7.13 0.8.2 (2021-08-27)

- **New functionality in `dev.labjack.LabJack`:**
  - configure clock and send timed pulse sequences
  - set DAC/analog output voltage
- Bugfix: ignore random bits sent by to `dev.newport.NewportSMC100PP` controller during start-up/powering-up.

## 7.14 0.8.1 (2021-08-13)

- Add Python version check (min version error; max version warning).
- Daily checks for upstream dependencies compatibility and devel environment improvements.

## 7.15 0.8.0 (2021-07-02)

- TCP communication protocol.
- Lauda PRO RP 245 E circulation thermostat device over TCP.
- Pico Technology PT-104 Platinum Resistance Data Logger device as a wrapper of the Python bindings for the PicoSDK.
- In `com.visa.VisaCommunication`: periodic status polling when VISA/TCP keep alive connection is not supported by a host.

## 7.16 0.7.1 (2021-06-04)

- New `utils.validation` submodule with `validate_bool` and `validate_number` utilities extracted from internal use within a `dev.tiepie` subpackage.
- In `comm.serial.SerialCommunication`:
  - strict encoding errors handling strategy for subclasses,
  - user warning for a low communication timeout value.

## 7.17 0.7.0 (2021-05-25)

- The `dev.tiepie` module was splitted into a subpackage with, in particular, submodules for each of the device types – `oscilloscope`, `generator`, and `i2c` – and with backward-incompatible direct imports from the submodules.
- In `dev.technix`:
  - fixed communication crash on nested status byte query;
  - added enums for GET and SET register commands.
- Further minor logging improvements: added missing module level logger and removed some error logs in `except` blocks used for a flow control.
- In `examples/` folder renamed consistently all the examples.
- In API documentation: fix incorrect links mapping on inheritance diagrams.

## 7.18 0.6.1 (2021-05-08)

- In `dev.tiepie`:
  - dynamically set oscilloscope's channel limits in `OscilloscopeChannelParameterLimits`: `input_range` and `trigger_level_abs`, incl. update of latter on each change of `input_range` value of a `TiePieOscilloscopeChannelConfig` instances;
  - quick fix for opening of combined instruments by disabling `OscilloscopeParameterLimits`. `trigger_delay` (an advanced feature);
  - enable automatic devices detection to be able to find network devices with `TiePieOscilloscope`. `list_devices()`.
- Fix `examples/example_labjack.py`.
- Improved logging: consistently use module level loggers, and always log exception tracebacks.
- Improve API documentation: separate pages per modules, each with an inheritance diagram as an overview.

## 7.19 0.6.0 (2021-04-23)

- Technix capacitor charger using either serial connection or Telnet protocol.
- **Extensions, improvements and fixes in existing devices:**
  - In `dev.tiepie.TiePieOscilloscope`:
    - \* redesigned measurement start and data collection API, incl. time out argument, with no/infinite time out option;
    - \* trigger allows now a no/infinite time out;
    - \* record length and trigger level were fixed to accept, respectively, floating point and integer numbers;
    - \* fixed resolution validation bug;
  - `dev.heinzinger.HeinzingerDI` and `dev.rs_rto1024.RTO1024` instances are now resilient to multiple `stop()` calls.
  - In `dev.crylas.CryLasLaser`: default configuration timeout and polling period were adjusted;
  - Fixed PSI9080 example script.
- **Package and source code improvements:**
  - Update to backward-incompatible `pyvisa-py>=0.5.2`. Developers, do update your local development environments!
  - External libraries, like LibTiePie SDK or LJM Library, are now not installed by default; they are now extra installation options.
  - Added Python 3.9 support.
  - Improved number formatting in logs.
  - Typing improvements and fixes for `mypy>=0.800`.

## 7.20 0.5.0 (2020-11-11)

- TiePie USB oscilloscope, generator and I2C host devices, as a wrapper of the Python bindings for the LibTiePie SDK.
- a FuG Elektronik Power Supply (e.g. Capacitor Charger HCK) using the built-in ADDAT controller with the Probus V protocol over a serial connection
- All devices polling status or measurements use now a `dev.utils.Poller` utility class.
- **Extensions and improvements in existing devices:**
  - In `dev.rs_rto1024.RT01024`: added Channel state, scale, range, position and offset accessors, and measurements activation and read methods.
  - In `dev.sst_luminox.Luminox`: added querying for all measurements in polling mode, and made output mode activation more robust.
  - In `dev.newport.NewportSMC100PP`: an error-prone `wait_until_move_finished` method of replaced by a fixed waiting time, device operations are now robust to a power supply cut, and device restart is not required to apply a start configuration.
- **Other minor improvements:**
  - Single failure-safe starting and stopping of devices sequenced via `dev.base.DeviceSequenceMixin`.
  - Moved `read_text_nonempty` up to `comm.serial.SerialCommunication`.
  - Added development Dockerfile.
  - Updated package and development dependencies: `pymodbus`, `pytest-mock`.

## 7.21 0.4.0 (2020-07-16)

- **Significantly improved new Supercube device controller:**
  - more robust error-handling,
  - status polling with generic Poller helper,
  - messages and status boards.
  - tested with a physical device,
- Improved OPC UA client wrapper, with better error handling, incl. re-tries on `concurrent.futures.TimeoutError`.
- SST Luminox Oxygen sensor device controller.
- **Backward-incompatible changes:**
  - `CommunicationProtocol.access_lock` has changed type from `threading.Lock` to `threading.RLock`.
  - `ILS2T.relative_step` and `ILS2T.absolute_position` are now called, respectively, `ILS2T.write_relative_step` and `ILS2T.write_absolute_position`.
- **Minor bugfixes and improvements:**
  - fix use of max resolution in `Labjack.set_ain_resolution()`,
  - resolve ILS2T devices relative and absolute position setters race condition,

- added acoustic horn function in the 2015 Supercube.

- **Toolchain changes:**

- add Python 3.8 support,
- drop pytest-runner support,
- ensure compatibility with `labjack_ljm` 2019 version library.

## 7.22 0.3.5 (2020-02-18)

- Fix issue with reading integers from LabJack LJM Library (device's product ID, serial number etc.)
- Fix development requirements specification (tox version).

## 7.23 0.3.4 (2019-12-20)

- **New devices using serial connection:**

- Heinzinger Digital Interface I/II and a Heinzinger PNC power supply
- Q-switched Pulsed Laser and a laser attenuator from CryLas
- Newport SMC100PP single axis motion controller for 2-phase stepper motors
- Pfeiffer TPG controller (TPG 25x, TPG 26x and TPG 36x) for Compact pressure Gauges

- PEP 561 compatibility and related corrections for static type checking (now in CI)

- **Refactorings:**

- Protected non-thread safe read and write in communication protocols
- Device sequence mixin: start/stop, add/rm and lookup
- `.format()` to f-strings
- more enumerations and a quite some improvements of existing code

- Improved error docstrings (`:raises:` annotations) and extended tests for errors.

## 7.24 0.3.3 (2019-05-08)

- Use PyPI labjack-ljm (no external dependencies)

## 7.25 0.3.2 (2019-05-08)

- INSTALLATION.rst with LJMPython prerequisite info

## 7.26 0.3.1 (2019-05-02)

- readthedocs.org support

## 7.27 0.3 (2019-05-02)

- Prevent an automatic close of VISA connection when not used.
- Rhode & Schwarz RTO 1024 oscilloscope using VISA interface over `TCP::INSTR`.
- Extended tests incl. messages sent to devices.
- Added Supercube device using an OPC UA client
- Added Supercube 2015 device using an OPC UA client (for interfacing with old system version)

## 7.28 0.2.1 (2019-04-01)

- Fix issue with LJMPython not being installed automatically with setuptools.

## 7.29 0.2.0 (2019-03-31)

- LabJack LJM Library communication wrapper and LabJack device.
- Modbus TCP communication protocol.
- Schneider Electric ILS2T stepper motor drive device.
- Elektro-Automatik PSI9000 current source device and VISA communication wrapper.
- Separate configuration classes for communication protocols and devices.
- Simple experiment manager class.

## 7.30 0.1.0 (2019-02-06)

- Communication protocol base and serial communication implementation.
- Device base and MBW973 implementation.



---

**CHAPTER  
EIGHT**

---

**INDICES AND TABLES**

- genindex
- modindex
- search



## PYTHON MODULE INDEX

### h

hvl\_ccb, 185  
hvl\_ccb.comm, 35  
hvl\_ccb.comm.base, 11  
hvl\_ccb.comm.labjack\_ljm, 16  
hvl\_ccb.comm.modbus\_tcp, 19  
hvl\_ccb.comm.opc, 21  
hvl\_ccb.comm.serial, 24  
hvl\_ccb.comm.tcp, 28  
hvl\_ccb.comm.telnet, 30  
hvl\_ccb.comm.visa, 32  
hvl\_ccb.configuration, 181  
hvl\_ccb.dev, 174  
hvl\_ccb.dev.base, 72  
hvl\_ccb.dev.crylas, 75  
hvl\_ccb.dev.cube, 48  
hvl\_ccb.dev.cube.base, 36  
hvl\_ccb.dev.cube.constants, 41  
hvl\_ccb.dev.cube.picube, 45  
hvl\_ccb.dev.ea\_psi9000, 85  
hvl\_ccb.dev.fluke884x, 57  
hvl\_ccb.dev.fluke884x.base, 48  
hvl\_ccb.dev.fluke884x.constants, 53  
hvl\_ccb.dev.fluke884x.ranges, 54  
hvl\_ccb.dev.fug, 90  
hvl\_ccb.dev.heinzinger, 103  
hvl\_ccb.dev.highland\_t560, 62  
hvl\_ccb.dev.highland\_t560.base, 57  
hvl\_ccb.dev.highland\_t560.channel, 59  
hvl\_ccb.dev.highland\_t560.device, 60  
hvl\_ccb.dev.labjack, 110  
hvl\_ccb.dev.lauda, 118  
hvl\_ccb.dev.mbw973, 125  
hvl\_ccb.dev.newport, 129  
hvl\_ccb.dev.pfeiffer\_tpg, 144  
hvl\_ccb.dev.rs\_rto1024, 150  
hvl\_ccb.dev.se\_ilis2t, 159  
hvl\_ccb.dev.sst\_luminox, 166  
hvl\_ccb.dev.technix, 68  
hvl\_ccb.dev.technix.base, 63  
hvl\_ccb.dev.technix.device, 66  
hvl\_ccb.dev.utils, 171

hvl\_ccb.dev.visa, 172  
hvl\_ccb.error, 183  
hvl\_ccb.experiment\_manager, 183  
hvl\_ccb.utils, 181  
hvl\_ccb.utils.conversion, 178  
hvl\_ccb.utils.conversion.map\_range, 174  
hvl\_ccb.utils.conversion.sensor, 175  
hvl\_ccb.utils.conversion.unit, 176  
hvl\_ccb.utils.conversion.utils, 177  
hvl\_ccb.utils.enum, 178  
hvl\_ccb.utils.typing, 180  
hvl\_ccb.utils.validation, 180



# INDEX

## A

A (*HeinzingerPNC.UnitCurrent attribute*), 107  
ABSOLUTE\_POSITION (*ILS2T.ActionsPtp attribute*), 159  
AC (*NewportConfigCommands attribute*), 129  
AC\_100KV (*PowerSetup attribute*), 44  
AC\_150KV (*PowerSetup attribute*), 44  
AC\_200KV (*PowerSetup attribute*), 44  
AC\_50KV (*PowerSetup attribute*), 44  
ac\_current\_range (*Fluke8845a attribute*), 49  
ac\_voltage\_range (*Fluke8845a attribute*), 49  
acceleration (*NewportSMC100PPConfig attribute*), 138  
ACCESS\_ENABLE (*ILS2TRegAddr attribute*), 165  
access\_lock (*CommunicationProtocol attribute*), 14  
ACCurrentRange (class in *hvl\_ccb.dev.fluke884x.ranges*), 54  
ACTION\_JOG\_VALUE (*ILS2T attribute*), 159  
activate\_clock\_output() (*T560 method*), 60  
activate\_measurements() (*RTO1024 method*), 151  
activate\_output() (*Luminox method*), 167  
activate\_remote\_mode() (*Fluke8845a method*), 49  
ACTIVE (*CryLasLaser.AnswersStatus attribute*), 78  
active\_alarms() (*BaseCube method*), 36  
ACTIVE\_HIGH (*Polarity attribute*), 58  
ACTIVE\_LOW (*Polarity attribute*), 58  
actualsetvalue (*FuGProbusVSetRegisters property*), 100  
ACVoltageRange (class in *hvl\_ccb.dev.fluke884x.ranges*), 55  
adc\_mode (*FuGProbusVMonitorRegisters property*), 99  
add\_device() (*DeviceSequenceMixin method*), 73  
add\_device() (*ExperimentManager method*), 184  
ADDR\_INCORRECT (*NewportSMC100PPSerialCommunication.ControllerBank attribute*), 139  
address (*NewportSMC100PPConfig attribute*), 138  
address (*VisaCommunicationConfig property*), 34  
address() (*VisaCommunicationConfig.InterfaceType method*), 33  
ADMODE (*FuGProbusIVCommands attribute*), 96  
all\_measurements (*LuminoxMeasurementType attribute*), 168

all\_measurements\_types() (*LuminoxMeasurementType class method*), 169  
ANALOG (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 122  
analog\_control (*FuGProbusVDIRegisters property*), 98  
ANY (*LabJack.DeviceType attribute*), 112  
ANY (*LJMCommunicationConfig.ConnectionType attribute*), 17  
ANY (*LJMCommunicationConfig.DeviceType attribute*), 17  
ApertureRange (class in *hvl\_ccb.dev.fluke884x.ranges*), 55  
AsyncCommunicationProtocol (class in *hvl\_ccb.comm.base*), 11  
AsyncCommunicationProtocolConfig (class in *hvl\_ccb.comm.base*), 13  
ATM (*Pressure attribute*), 176  
ATMOSPHERE (*Pressure attribute*), 176  
attenuation (*CryLasAttenuator property*), 75  
AUTO (*EarthingStickOperatingStatus attribute*), 42  
AUTO (*LaudaProRp245eConfig.OperationModeEnum attribute*), 122  
AUTO (*RTO1024.TriggerModes attribute*), 151  
auto\_install\_mode (*T560 property*), 60  
auto\_install\_mode (*T560Config attribute*), 61  
auto\_laser\_on (*CryLasLaserConfig attribute*), 82  
AutoInstallMode (class in *hvl\_ccb.dev.highland\_t560.base*), 57  
AutoNumberNameEnum (class in *hvl\_ccb.utils.enum*), 178

## B

BA (*NewportConfigCommands attribute*), 129  
backlash\_compensation (*NewportSMC100PPConfig attribute*), 138  
backup\_waveform() (*RTO1024 method*), 151  
BAR (*Pressure attribute*), 176  
barometric\_pressure (*LuminoxMeasurementType attribute*), 169  
base\_velocity (*NewportSMC100PPConfig attribute*), 138  
BaseCube (class in *hvl\_ccb.dev.cube.base*), 36

BaseCubeConfiguration (class *in* C  
    *hvl\_ccb.dev(cube.base)*, 39  
BaseCubeOpcUaCommunication (class *in* C  
    *hvl\_ccb.dev(cube.base)*, 40  
BaseCubeOpcUaCommunicationConfig (class *in* C  
    *hvl\_ccb.dev(cube.base)*, 40  
BATH\_TEMP (*LaudaProRp245eCommand attribute*), 120  
baudrate (*CryLasAttenuatorSerialCommunicationConfig attribute*), 77  
baudrate (*CryLasLaserSerialCommunicationConfig attribute*), 84  
baudrate (*FuGSerialCommunicationConfig attribute*), 102  
baudrate (*HeinzingerSerialCommunicationConfig attribute*), 109  
baudrate (*LuminoxSerialCommunicationConfig attribute*), 170  
baudrate (*MBW973SerialCommunicationConfig attribute*), 128  
baudrate (*NewportSMC100PPSerialCommunicationConfig attribute*), 142  
baudrate (*PfeifferTPGSerialCommunicationConfig attribute*), 149  
baudrate (*SerialCommunicationConfig attribute*), 26  
BH (*NewportConfigCommands attribute*), 129  
board (*VisaCommunicationConfig attribute*), 34  
BoolEnum (class *in hvl\_ccb.utils.enum*), 178  
breakdown\_detection\_active (*BaseCube property*), 36  
breakdown\_detection\_reset() (*BaseCube method*), 36  
breakdown\_detection\_triggered (*BaseCube property*), 36  
bufsize (*TcpCommunicationConfig attribute*), 29  
build\_str() (*LaudaProRp245eCommand method*), 121  
BUS (*TriggerSource attribute*), 54  
bytesize (*CryLasAttenuatorSerialCommunicationConfig attribute*), 77  
bytesize (*CryLasLaserSerialCommunicationConfig attribute*), 84  
bytesize (*FuGSerialCommunicationConfig attribute*), 102  
bytesize (*HeinzingerSerialCommunicationConfig attribute*), 109  
bytesize (*LuminoxSerialCommunicationConfig attribute*), 170  
bytesize (*MBW973SerialCommunicationConfig attribute*), 128  
bytesize (*NewportSMC100PPSerialCommunicationConfig attribute*), 142  
bytesize (*PfeifferTPGSerialCommunicationConfig attribute*), 149  
Bytesize (*SerialCommunicationConfig attribute*), 26  
bytesize (*SerialCommunicationConfig attribute*), 26  
    *in* C  
        *C (LabJack.TemperatureUnit attribute)*, 113  
        *C (LabJack.ThermocoupleType attribute)*, 113  
        *C (Temperature attribute)*, 177  
        *calibration\_factor (CryLasLaserConfig attribute)*, 82  
        *calibration\_factor (LEM4000S attribute)*, 175  
        *calibration\_mode (FuGProbusVDIRegisters property)*, 98  
        *cc\_mode (FuGProbusVDIRegisters property)*, 98  
        CCBErro, 183  
        cee16\_socket (*BaseCube property*), 36  
        CELSIUS (*Temperature attribute*), 177  
        ch\_a (*T560 property*), 60  
        ch\_b (*T560 property*), 60  
        ch\_c (*T560 property*), 60  
        ch\_d (*T560 property*), 60  
        check\_for\_error() (*NewportSMC100PPSerialCommunication method*), 140  
        check\_generic\_type() (*in module hvl\_ccb.utils.typing*), 180  
        check\_master\_slave\_config() (*PSI9000 method*), 86  
        chunk\_size (*VisaCommunicationConfig attribute*), 34  
        clean\_values() (*AsyncCommunicationProtocolConfig method*), 13  
        clean\_values() (*BaseCubeConfiguration method*), 39  
        clean\_values() (*CryLasAttenuatorConfig method*), 76  
        clean\_values() (*CryLasLaserConfig method*), 82  
        clean\_values() (*EmptyConfig method*), 74, 182  
        clean\_values() (*Fluke8845aConfig method*), 51  
        clean\_values() (*FuGConfig method*), 92  
        clean\_values() (*HeinzingerConfig method*), 104  
        clean\_values() (*ILS2TConfig method*), 163  
        clean\_values() (*LaudaProRp245eConfig method*), 122  
        clean\_values() (*LaudaProRp245eTcpCommunicationConfig method*), 124  
        clean\_values() (*LJMCommunicationConfig method*), 18  
        clean\_values() (*LuminoxConfig method*), 168  
        clean\_values() (*MBW973Config method*), 127  
        clean\_values() (*ModbusTcpCommunicationConfig method*), 20  
        clean\_values() (*NewportSMC100PPConfig method*), 138  
        clean\_values() (*OpcUaCommunicationConfig method*), 22  
        clean\_values() (*PfeifferTPGConfig method*), 147  
        clean\_values() (*PICubeConfiguration method*), 46  
        clean\_values() (*PSI9000Config method*), 88

**clean\_values()** (*SerialCommunicationConfig method*), 26  
**clean\_values()** (*T560Config method*), 61  
**clean\_values()** (*TcpCommunicationConfig method*), 29  
**clean\_values()** (*TechnixConfig method*), 67  
**clean\_values()** (*TelnetCommunicationConfig method*), 31  
**clean\_values()** (*VisaCommunicationConfig method*), 34  
**clear\_display\_message()** (*Fluke8845a method*), 49  
**clear\_error\_queue()** (*Fluke8845a method*), 49  
**Client** (*class in hvl\_ccb.comm.opc*), 21  
**CLOSE** (*EarthingStickOperation attribute*), 42  
**close()** (*CommunicationProtocol method*), 14  
**close()** (*LaudaProRp245eTcpCommunication method*), 123  
**close()** (*LJMCommunication method*), 16  
**close()** (*ModbusTcpCommunication method*), 19  
**close()** (*NullCommunicationProtocol method*), 14  
**close()** (*OpcUaCommunication method*), 21  
**close()** (*SerialCommunication method*), 24  
**close()** (*Tcp method*), 28  
**close()** (*TelnetCommunication method*), 30  
**close()** (*VisaCommunication method*), 32  
**close\_shutter()** (*CryLasLaser method*), 79  
**CLOSED** (*CryLasLaser.AnswersShutter attribute*), 78  
**CLOSED** (*CryLasLaserShutterStatus attribute*), 85  
**CLOSED** (*DoorStatus attribute*), 42  
**CLOSED** (*EarthingStickStatus attribute*), 43  
**CMD\_EXEC\_ERROR** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_CC** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_CONFIGURATION** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_DISABLE** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_HOMING** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_MOVING** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_NOT\_REFERENCED** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140

**CMD\_NOT\_ALLOWED\_PP** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMD\_NOT\_ALLOWED\_READY** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**CMR** (*PfeifferTPG.SensorTypes attribute*), 145  
**CODE\_OR\_ADDR\_INVALID** (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**com** (*SingleCommDevice property*), 74  
**COM\_TIME\_OUT** (*LaudaProRp245eCommand attribute*), 120  
**com\_time\_out** (*LaudaProRp245eConfig attribute*), 122  
**COM\_TIMEOUT** (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**command** (*LuminoxMeasurementType property*), 169  
**COMMAND** (*TriggerMode attribute*), 59  
**command()** (*FuGProbusIV method*), 95  
**communication\_channel** (*TechnixConfig attribute*), 67  
**CommunicationError**, 14  
**CommunicationProtocol** (*class in hvl\_ccb.comm.base*), 14  
**config** (*ConfigurationMixin property*), 181  
**CONFIG** (*FuGProbusVRegisterGroups attribute*), 99  
**CONFIG** (*NewportSMC100PP.StateMessages attribute*), 131  
**CONFIG** (*NewportStates attribute*), 143  
**config\_cls()** (*AsyncCommunicationProtocol static method*), 11  
**config\_cls()** (*BaseCube static method*), 37  
**config\_cls()** (*BaseCubeOpcUaCommunication static method*), 40  
**config\_cls()** (*ConfigurationMixin static method*), 181  
**config\_cls()** (*CryLasAttenuator static method*), 75  
**config\_cls()** (*CryLasAttenuatorSerialCommunication static method*), 77  
**config\_cls()** (*CryLasLaser static method*), 79  
**config\_cls()** (*CryLasLaserSerialCommunication static method*), 83  
**config\_cls()** (*Device static method*), 72  
**config\_cls()** (*Fluke8845a static method*), 49  
**config\_cls()** (*Fluke8845aTelnetCommunication static method*), 51  
**config\_cls()** (*FuGProbusIV static method*), 96  
**config\_cls()** (*FuGSerialCommunication static method*), 101  
**config\_cls()** (*HeinzingerDI static method*), 105  
**config\_cls()** (*HeinzingerSerialCommunication static method*), 108  
**config\_cls()** (*ILS2T static method*), 160  
**config\_cls()** (*ILS2TModbusTcpCommunication static method*), 164  
**config\_cls()** (*LaudaProRp245e static method*), 118

**config\_cls()** (*LaudaProRp245eTcpCommunication static method*), 123  
**config\_cls()** (*LJMCommunication static method*), 16  
**config\_cls()** (*Luminox static method*), 167  
**config\_cls()** (*LuminoxSerialCommunication static method*), 170  
**config\_cls()** (*MBW973 static method*), 125  
**config\_cls()** (*MBW973SerialCommunication static method*), 128  
**config\_cls()** (*ModbusTcpCommunication static method*), 19  
**config\_cls()** (*NewportSMC100PP static method*), 132  
**config\_cls()** (*NewportSMC100PPSerialCommunication static method*), 140  
**config\_cls()** (*NullCommunicationProtocol static method*), 15  
**config\_cls()** (*OpcUaCommunication static method*), 21  
**config\_cls()** (*PfeifferTPG static method*), 145  
**config\_cls()** (*PfeifferTPGSerialCommunication static method*), 148  
**config\_cls()** (*PICube static method*), 45  
**config\_cls()** (*PICubeOpcUaCommunication static method*), 47  
**config\_cls()** (*PSI9000 static method*), 86  
**config\_cls()** (*PSI9000VisaCommunication static method*), 89  
**config\_cls()** (*RTO1024 static method*), 151  
**config\_cls()** (*RTO1024VisaCommunication static method*), 158  
**config\_cls()** (*SerialCommunication static method*), 25  
**config\_cls()** (*SyncCommunicationProtocol static method*), 15  
**config\_cls()** (*T560 static method*), 60  
**config\_cls()** (*T560Communication static method*), 58  
**config\_cls()** (*Tcp static method*), 28  
**config\_cls()** (*Technix static method*), 66  
**config\_cls()** (*TechnixSerialCommunication static method*), 63  
**config\_cls()** (*TechnixTelnetCommunication static method*), 64  
**config\_cls()** (*TelnetCommunication static method*), 30  
**config\_cls()** (*VisaCommunication static method*), 32  
**config\_cls()** (*VisaDevice static method*), 172  
**config\_high\_pulse()** (*LabJack method*), 113  
**config\_status** (*FuG property*), 90  
**configdataclass()** (*in module hvl\_ccb.configuration*), 182  
**configuration\_save\_json()** (*ConfigurationMixin method*), 181  
**ConfigurationMixin** (*class in hvl\_ccb.configuration*), 181  
**ConfigurationValueWarning**, 182  
**connection\_type** (*LJMCommunicationConfig attribute*), 18  
**CONT\_MODE** (*LaudaProRp245eCommand attribute*), 120  
**continue\_ramp()** (*LaudaProRp245e method*), 118  
**control\_mode** (*LaudaProRp245eConfig attribute*), 122  
**CONVERSION** (*LEM4000S attribute*), 175  
**convert()** (*LEM4000S method*), 175  
**convert()** (*LMT70A method*), 176  
**convert()** (*Sensor method*), 176  
**convert()** (*Unit class method*), 177  
**convert\_to\_range1()** (*MapRanges method*), 175  
**convert\_to\_range2()** (*MapRanges method*), 175  
**convert\_value\_to\_str()** (*in module hvl\_ccb.utils.conversion.utils*), 177  
**ConvertibleTypes** (*in module hvl\_ccb.utils.typing*), 180  
**COOLOFF** (*LaudaProRp245eConfig.OperationModeEnum attribute*), 122  
**COOLON** (*LaudaProRp245eConfig.OperationModeEnum attribute*), 122  
**CR** (*FuGTerminators attribute*), 103  
**create\_serial\_port()** (*SerialCommunicationConfig method*), 26  
**create\_telnet()** (*TelnetCommunicationConfig method*), 31  
**CRLF** (*FuGTerminators attribute*), 103  
**CryLasAttenuator** (*class in hvl\_ccb.dev.crylas*), 75  
**CryLasAttenuatorConfig** (*class in hvl\_ccb.dev.crylas*), 76  
**CryLasAttenuatorError**, 77  
**CryLasAttenuatorSerialCommunication** (*class in hvl\_ccb.dev.crylas*), 77  
**CryLasAttenuatorSerialCommunicationConfig** (*class in hvl\_ccb.dev.crylas*), 77  
**CryLasLaser** (*class in hvl\_ccb.dev.crylas*), 78  
**CryLasLaser.AnswersShutter** (*class in hvl\_ccb.dev.crylas*), 78  
**CryLasLaser.AnswersStatus** (*class in hvl\_ccb.dev.crylas*), 78  
**CryLasLaser.LaserStatus** (*class in hvl\_ccb.dev.crylas*), 79  
**CryLasLaser.RepetitionRates** (*class in hvl\_ccb.dev.crylas*), 79  
**CryLasLaserConfig** (*class in hvl\_ccb.dev.crylas*), 81  
**CryLasLaserError**, 82  
**CryLasLaserNotReadyError**, 82  
**CryLasLaserPoller** (*class in hvl\_ccb.dev.crylas*), 83  
**CryLasLaserSerialCommunication** (*class in hvl\_ccb.dev.crylas*), 83  
**CryLasLaserSerialCommunicationConfig** (*class in hvl\_ccb.dev.crylas*), 84  
**CryLasLaserShutterStatus** (*class in hvl\_ccb.dev.crylas*), 85  
**CubeEarthingStickOperationError**, 41

**CubeError**, 41  
**CubeRemoteControlError**, 42  
**CubeStatusChangeError**, 42  
**CubeStopError**, 42  
**current (FuG property)**, 91  
**CURRENT (FuGProbusIVCommands attribute)**, 96  
**CURRENT (FuGReadbackChannels attribute)**, 101  
**current (Technix property)**, 66  
**CURRENT\_AC (MeasurementFunction attribute)**, 53  
**CURRENT\_DC (MeasurementFunction attribute)**, 53  
**current\_filter (Fluke8845a attribute)**, 49  
**current\_lower\_limit (PSI9000Config attribute)**, 88  
**current\_monitor (FuG property)**, 91  
**current\_primary (PICube property)**, 45  
**current\_upper\_limit (PSI9000Config attribute)**, 88  
**CurrentPosition** (New-  
portSMC100PPConfig.HomeSearch attribute), 138  
**cv\_mode (FuGProbusVDIRegisters property)**, 98

**D**

**datachange\_notification()** (*OpcUaSubHandler method*), 24  
**date\_of\_manufacture (LuminoxMeasurementType attribute)**, 169  
**datetime\_to\_opc () (BaseCube class method)**, 37  
**DC\_140KV (PowerSetup attribute)**, 44  
**DC\_280KV (PowerSetup attribute)**, 44  
**dc\_current\_range (Fluke8845a attribute)**, 49  
**dc\_voltage\_range (Fluke8845a attribute)**, 49  
**DC\_VOLTAGE\_TOO\_LOW** (New-  
portSMC100PP.MotorErrors attribute), 130  
**DCCurrentRange** (class in hvl\_ccb.dev.fluke884x.ranges), 55  
**DCVoltageRange** (class in hvl\_ccb.dev.fluke884x.ranges), 55  
**default\_com\_cls () (BaseCube static method)**, 37  
**default\_com\_cls () (CryLasAttenuator static method)**, 75  
**default\_com\_cls () (CryLasLaser static method)**, 79  
**default\_com\_cls () (Fluke8845a static method)**, 49  
**default\_com\_cls () (FuGProbusIV static method)**, 96  
**default\_com\_cls () (HeinzingerDI static method)**, 105  
**default\_com\_cls () (ILS2T static method)**, 160  
**default\_com\_cls () (LabJack static method)**, 113  
**default\_com\_cls () (LaudaProRp245e static method)**, 118  
**default\_com\_cls () (Luminox static method)**, 167  
**default\_com\_cls () (MBW973 static method)**, 125  
**default\_com\_cls () (NewportSMC100PP static method)**, 132  
**default\_com\_cls () (PfeifferTPG static method)**, 145  
**default\_com\_cls () (PICube static method)**, 45

**default\_com\_cls () (PSI9000 static method)**, 86  
**default\_com\_cls () (RTO1024 static method)**, 151  
**default\_com\_cls () (SingleCommDevice static method)**, 74  
**default\_com\_cls () (T560 static method)**, 61  
**default\_com\_cls () (Technix method)**, 66  
**default\_com\_cls () (VisaDevice static method)**, 172  
**DEFAULT\_IO\_SCANNING\_CONTROL\_VALUES (ILS2T attribute)**, 159  
**default\_n\_attempts\_read\_text\_nonempty (Async-CommunicationProtocolConfig attribute)**, 13  
**default\_n\_attempts\_read\_text\_nonempty (FuGSerialCommunicationConfig attribute)**, 102  
**default\_n\_attempts\_read\_text\_nonempty (HeinzingerSerialCommunicationConfig attribute)**, 109  
**default\_number\_of\_recordings (HeinzingerConfig attribute)**, 104  
**Device** (class in hvl\_ccb.dev.base), 72  
**DEVICE\_TYPE (LaudaProRp245eCommand attribute)**, 121  
**device\_type (LJMCommunicationConfig attribute)**, 18  
**DeviceError**, 72  
**DeviceExistingError**, 72  
**DeviceFailuresError**, 72  
**devices\_failed\_start (DeviceSequenceMixin attribute)**, 73  
**devices\_failed\_start (ExperimentManager attribute)**, 184  
**devices\_failed\_stop (DeviceSequenceMixin attribute)**, 73  
**devices\_failed\_stop (ExperimentManager attribute)**, 184  
**DeviceSequenceMixin** (class in hvl\_ccb.dev.base), 72  
**di (FuG property)**, 91  
**digital\_control (FuGProbusVDIRegisters property)**, 98  
**DIOChannel (LabJack attribute)**, 112  
**DIODE (MeasurementFunction attribute)**, 53  
**DISABLE (NewportStates attribute)**, 143  
**disable () (ILS2T method)**, 160  
**DISABLE\_FROM\_JOGGING** (New-  
portSMC100PP.StateMessages attribute), 131  
**DISABLE\_FROM\_MOVING** (New-  
portSMC100PP.StateMessages attribute), 131  
**DISABLE\_FROM\_READY** (New-  
portSMC100PP.StateMessages attribute), 131  
**disable\_pulses () (LabJack method)**, 114  
**DisableEspStageCheck** (New-  
portSMC100PPConfig.EspStageConfig attribute), 137

disarm_trigger() ( <i>T560 method</i> ), 61			
disconnect() ( <i>Client method</i> ), 21			
DISPLACEMENT_OUT_OF_LIMIT	(New-	EarthingStickOperation	(class
<i>portSMC100PPSerialCommunication.ControllerErrors</i>		<i>hvl_ccb.dev(cube.constants)</i> ), 42	in
attribute), 140			
display_enable ( <i>Fluke8845a property</i> ), 49		EarthingStickStatus	(class
DISPLAY_MAX_LENGTH ( <i>Fluke8845a attribute</i> ), 49		<i>hvl_ccb.dev(cube.constants)</i> ), 43	in
display_message ( <i>Fluke8845a property</i> ), 49		EEPROM_ACCESS_ERROR	(New-
display_message_board() ( <i>BaseCube method</i> ), 37		<i>portSMC100PPSerialCommunication.ControllerErrors</i>	
display_status_board() ( <i>BaseCube method</i> ), 37		attribute), 140	
do_iосcanning_write() ( <i>ILS2T method</i> ), 160		EIGHT	( <i>HeinzingerConfig.RecordingsEnum attribute</i> ),
door_1_status ( <i>BaseCube attribute</i> ), 37		104	
door_2_status ( <i>BaseCube attribute</i> ), 37		EIGHTBITS	( <i>SerialCommunicationBytesize attribute</i> ), 25
door_3_status ( <i>BaseCube attribute</i> ), 37		EmptyConfig	( <i>class in hvl_ccb.configuration</i> ), 182
DoorStatus ( <i>class in hvl_ccb.dev(cube.constants)</i> ), 42		EmptyConfig	( <i>class in hvl_ccb.dev.base</i> ), 74
<b>E</b>		enable()	( <i>ILS2T method</i> ), 161
E ( <i>LabJack.ThermocoupleType attribute</i> ), 113		enable_clock()	( <i>LabJack method</i> ), 114
E0 ( <i>FuGErrorcodes attribute</i> ), 93		EnableEspStageCheck	(New-
E1 ( <i>FuGErrorcodes attribute</i> ), 93		<i>portSMC100PPConfig.EspStageConfig attribute</i> ), 138	
E10 ( <i>FuGErrorcodes attribute</i> ), 93		encoding	( <i>AsyncCommunicationProtocolConfig attribute</i> ), 13
E100 ( <i>FuGErrorcodes attribute</i> ), 93		encoding	( <i>NewportSMC100PPSerialCommunicationConfig attribute</i> ), 142
E106 ( <i>FuGErrorcodes attribute</i> ), 94		encoding_error_handling	( <i>AsyncCommunicationProtocolConfig attribute</i> ), 13
E11 ( <i>FuGErrorcodes attribute</i> ), 94		encoding_error_handling	( <i>NewportSMC100PPSerialCommunicationConfig attribute</i> ), 142
E115 ( <i>FuGErrorcodes attribute</i> ), 94		EndOfRunSwitch	(New-
E12 ( <i>FuGErrorcodes attribute</i> ), 94		<i>portSMC100PPConfig.HomeSearch attribute</i> ), 138	
E125 ( <i>FuGErrorcodes attribute</i> ), 94		EndOfRunSwitch_and_Index	(New-
E13 ( <i>FuGErrorcodes attribute</i> ), 94		<i>portSMC100PPConfig.HomeSearch attribute</i> ), 138	
E135 ( <i>FuGErrorcodes attribute</i> ), 94		endpoint_name	( <i>BaseCubeOpcUaCommunicationConfig attribute</i> ), 41
E14 ( <i>FuGErrorcodes attribute</i> ), 94		endpoint_name	( <i>OpcUaCommunicationConfig attribute</i> ), 23
E145 ( <i>FuGErrorcodes attribute</i> ), 94		endpoint_name	( <i>PICubeOpcUaCommunicationConfig attribute</i> ), 47
E15 ( <i>FuGErrorcodes attribute</i> ), 94		ERROR	( <i>DoorStatus attribute</i> ), 42
E155 ( <i>FuGErrorcodes attribute</i> ), 94		ERROR	( <i>EarthingStickStatus attribute</i> ), 43
E16 ( <i>FuGErrorcodes attribute</i> ), 94		ERROR	( <i>ExperimentStatus attribute</i> ), 185
E165 ( <i>FuGErrorcodes attribute</i> ), 94		ERROR	( <i>SafetyStatus attribute</i> ), 44
E2 ( <i>FuGErrorcodes attribute</i> ), 94		errorcode	( <i>FuGError attribute</i> ), 93
E206 ( <i>FuGErrorcodes attribute</i> ), 94		ESP_STAGE_NAME_INVALID	(New-
E306 ( <i>FuGErrorcodes attribute</i> ), 94		<i>portSMC100PPSerialCommunication.ControllerErrors attribute</i> ), 140	
E4 ( <i>FuGErrorcodes attribute</i> ), 94		ETH	( <i>LaudaProRp245eConfig.ExtControlModeEnum attribute</i> ), 122
E5 ( <i>FuGErrorcodes attribute</i> ), 94		ETHERNET	( <i>LJMCommunicationConfig.ConnectionType attribute</i> ), 17
E504 ( <i>FuGErrorcodes attribute</i> ), 94		EVEN	( <i>SerialCommunicationParity attribute</i> ), 27
E505 ( <i>FuGErrorcodes attribute</i> ), 94		event_notification()	( <i>OpcUaSubHandler method</i> ), 24
E6 ( <i>FuGErrorcodes attribute</i> ), 94		EXECUTE	( <i>FuGProbusIVCommands attribute</i> ), 96
E666 ( <i>FuGErrorcodes attribute</i> ), 95			
E7 ( <i>FuGErrorcodes attribute</i> ), 95			
E8 ( <i>FuGErrorcodes attribute</i> ), 95			
E9 ( <i>FuGErrorcodes attribute</i> ), 95			
earthing_rod_1_status ( <i>BaseCube attribute</i> ), 37			
earthing_rod_2_status ( <i>BaseCube attribute</i> ), 37			
earthing_rod_3_status ( <i>BaseCube attribute</i> ), 37			
EarthingRodStatus	(class		
<i>hvl_ccb.dev(cube.constants)</i> ), 42	in		
EarthingStickOperatingStatus	(class		
<i>hvl_ccb.dev(cube.constants)</i> ), 42	in		

**execute\_absolute\_position()** (*ILS2T method*), 161  
**execute\_on\_x** (*FuGProbusVConfigRegisters property*), 97  
**execute\_relative\_step()** (*ILS2T method*), 161  
**EXECUTEONX** (*FuGProbusIVCommands attribute*), 96  
**exit\_configuration()** (*NewportSMC100PP method*), 132  
**exit\_configuration\_wait\_sec** (*NewportSMC100PPConfig attribute*), 138  
**EXPERIMENT\_BLOCKED** (*EarthingRodStatus attribute*), 42  
**EXPERIMENT\_READY** (*EarthingRodStatus attribute*), 42  
**ExperimentError**, 183  
**ExperimentManager** (class in *hvl\_ccb.experiment\_manager*), 183  
**ExperimentStatus** (class in *hvl\_ccb.experiment\_manager*), 185  
**EXPT100** (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 122  
**EXT\_FALLING\_EDGE** (*TriggerMode attribute*), 59  
**EXT\_RISING\_EDGE** (*TriggerMode attribute*), 59  
**EXTERNAL** (*TriggerSource attribute*), 54  
**EXTERNAL\_SOURCE** (*PowerSetup attribute*), 44  
**EXTERNAL\_TEMP** (*LaudaProRp245eCommand attribute*), 121

**F**

**F** (*LabJack.TemperatureUnit attribute*), 113  
**F** (*Temperature attribute*), 177  
**FARENHEIT** (*Temperature attribute*), 177  
**failures** (*DeviceFailuresError attribute*), 72  
**FAST** (*ILS2T.Ref16Jog attribute*), 159  
**FAST\_FILTER** (*FilterRange attribute*), 56  
**fetch()** (*Fluke8845a method*), 49  
**file\_copy()** (*RTO1024 method*), 151  
**FilterRange** (class in *hvl\_ccb.dev.fluke884x.ranges*), 56  
**finish()** (*ExperimentManager method*), 184  
**FINISHED** (*ExperimentStatus attribute*), 185  
**FINISHING** (*ExperimentStatus attribute*), 185  
**fire\_trigger()** (*T560 method*), 61  
**FIRMWARE** (*FuGReadbackChannels attribute*), 101  
**FIVE\_MHZ** (*LabJack.ClockFrequency attribute*), 112  
**FIVEBITS** (*SerialCommunicationBytesize attribute*), 25  
**FLT\_INFO** (*ILS2TRegAddr attribute*), 165  
**FLT\_MEM\_DEL** (*ILS2TRegAddr attribute*), 165  
**FLT\_MEM\_RESET** (*ILS2TRegAddr attribute*), 165  
**Fluke8845a** (class in *hvl\_ccb.dev.fluke884x.base*), 48  
**Fluke8845aCheckError**, 53  
**Fluke8845aConfig** (class in *hvl\_ccb.dev.fluke884x.base*), 51  
**Fluke8845aError**, 53  
**Fluke8845aTelnetCommunication** (class in *hvl\_ccb.dev.fluke884x.base*), 51  
**Fluke8845aTelnetCommunicationConfig** (class in *hvl\_ccb.dev.fluke884x.base*), 52  
**Fluke8845aUnknownCommandError**, 53  
**FOLLOWING\_ERROR** (*NewportSMC100PP.MotorErrors attribute*), 130  
**FOLLOWRAMP** (*FuGRampModes attribute*), 101  
**force\_value()** (*AsyncCommunicationProtocolConfig method*), 13  
**force\_value()** (*BaseCubeConfiguration method*), 39  
**force\_value()** (*BaseCubeOpcUaCommunicationConfig method*), 41  
**force\_value()** (*CryLasAttenuatorConfig method*), 76  
**force\_value()** (*CryLasAttenuatorSerialCommunicationConfig method*), 77  
**force\_value()** (*CryLasLaserConfig method*), 82  
**force\_value()** (*CryLasLaserSerialCommunicationConfig method*), 84  
**force\_value()** (*EmptyConfig method*), 74, 182  
**force\_value()** (*Fluke8845aConfig method*), 51  
**force\_value()** (*Fluke8845aTelnetCommunicationConfig method*), 52  
**force\_value()** (*FuGConfig method*), 92  
**force\_value()** (*FuGSerialCommunicationConfig method*), 102  
**force\_value()** (*HeinzingerConfig method*), 104  
**force\_value()** (*HeinzingerSerialCommunicationConfig method*), 109  
**force\_value()** (*ILS2TConfig method*), 163  
**force\_value()** (*ILS2TModbusTcpCommunicationConfig method*), 164  
**force\_value()** (*LaudaProRp245eConfig method*), 122  
**force\_value()** (*LaudaProRp245eTcpCommunicationConfig method*), 124  
**force\_value()** (*LJMCommunicationConfig method*), 18  
**force\_value()** (*LuminoxConfig method*), 168  
**force\_value()** (*LuminoxSerialCommunicationConfig method*), 170  
**force\_value()** (*MBW973Config method*), 127  
**force\_value()** (*MBW973SerialCommunicationConfig method*), 128  
**force\_value()** (*ModbusTcpCommunicationConfig method*), 20  
**force\_value()** (*NewportSMC100PPConfig method*), 138  
**force\_value()** (*NewportSMC100PPSerialCommunicationConfig method*), 142  
**force\_value()** (*OpcUaCommunicationConfig method*), 23  
**force\_value()** (*PfeifferTPGConfig method*), 147  
**force\_value()** (*PfeifferTPGSerialCommunicationConfig method*), 149  
**force\_value()** (*PICubeConfiguration method*), 46  
**force\_value()** (*PICubeOpcUaCommunicationConfig method*), 47

force_value()	( <i>PSI9000Config</i> method), 88		
force_value()	( <i>PSI9000VisaCommunicationConfig</i> method), 89		
force_value()	( <i>RTO1024Config</i> method), 157		
force_value()	( <i>RTO1024VisaCommunicationConfig</i> method), 158		
force_value()	( <i>SerialCommunicationConfig</i> method), 26		
force_value()	( <i>T560CommunicationConfig</i> method), 58		
force_value()	( <i>T560Config</i> method), 61		
force_value()	( <i>TcpCommunicationConfig</i> method), 29		
force_value()	( <i>TechnixConfig</i> method), 67		
force_value()	( <i>TechnixSerialCommunicationConfig</i> method), 64		
force_value()	( <i>TechnixTelnetCommunicationConfig</i> method), 65		
force_value()	( <i>TelnetCommunicationConfig</i> method), 31		
force_value()	( <i>VisaCommunicationConfig</i> method), 34		
force_value()	( <i>VisaDeviceConfig</i> method), 173		
FORTY_MHZ	( <i>LabJack.ClockFrequency</i> attribute), 112		
FOUR	( <i>HeinzingerConfig.RecordingsEnum</i> attribute), 104		
FOUR_HUNDRED_MILLI_AMPERE	( <i>ACCurrrentRange</i> attribute), 54		
FOUR_HUNDRED_MILLI_AMPERE	( <i>DCCurrentRange</i> attribute), 55		
FOUR_WIRE_RESISTANCE	( <i>MeasurementFunction</i> attribute), 53		
four_wire_resistance_range	( <i>Fluke8845a</i> attribute), 50		
FREERUN	( <i>RTO1024.TriggerModes</i> attribute), 151		
FREQUENCY	( <i>MeasurementFunction</i> attribute), 53		
frequency	( <i>PICube</i> property), 45		
frequency	( <i>T560</i> property), 61		
frequency_aperture	( <i>Fluke8845a</i> attribute), 50		
FRM	( <i>NewportConfigCommands</i> attribute), 129		
from_json()	( <i>ConfigurationMixin</i> class method), 182		
FRS	( <i>NewportConfigCommands</i> attribute), 129		
FuG	(class in <i>hvl_ccb.dev.fug</i> ), 90		
FuGConfig	(class in <i>hvl_ccb.dev.fug</i> ), 92		
FuGDigitalVal	(class in <i>hvl_ccb.dev.fug</i> ), 93		
FuGError	, 93		
FuGErrorcodes	(class in <i>hvl_ccb.dev.fug</i> ), 93		
FuGMonitorModes	(class in <i>hvl_ccb.dev.fug</i> ), 95		
FuGPolarities	(class in <i>hvl_ccb.dev.fug</i> ), 95		
FuGProbusIV	(class in <i>hvl_ccb.dev.fug</i> ), 95		
FuGProbusIVCommands	(class in <i>hvl_ccb.dev.fug</i> ), 96		
FuGProbusV	(class in <i>hvl_ccb.dev.fug</i> ), 97		
FuGProbusVConfigRegisters	(class in <i>hvl_ccb.dev.fug</i> ), 97		
FuGProbusVDIRegisters	(class in <i>hvl_ccb.dev.fug</i> ), 98		
FuGProbusVDORegisters	(class in <i>hvl_ccb.dev.fug</i> ), 99		
FuGProbusVMonitorRegisters	(class in <i>hvl_ccb.dev.fug</i> ), 99		
FuGProbusVRegisterGroups	(class in <i>hvl_ccb.dev.fug</i> ), 99		
FuGProbusVSetRegisters	(class in <i>hvl_ccb.dev.fug</i> ), 100		
FuGRampModes	(class in <i>hvl_ccb.dev.fug</i> ), 100		
FuReadbackChannels	(class in <i>hvl_ccb.dev.fug</i> ), 101		
FuGSerialCommunication	(class in <i>hvl_ccb.dev.fug</i> ), 101		
FuGSerialCommunicationConfig	(class in <i>hvl_ccb.dev.fug</i> ), 102		
FuGTerminators	(class in <i>hvl_ccb.dev.fug</i> ), 103		
<b>G</b>			
gate_mode	( <i>T560</i> property), 61		
gate_polarity	( <i>T560</i> property), 61		
GateMode	(class in <i>hvl_ccb.dev.hIGHLAND_t560.base</i> ), 57		
get_acceleration()	( <i>NewportSMC100PP</i> method), 132		
get_acquire_length()	( <i>RTO1024</i> method), 152		
get_ain()	( <i>LabJack</i> method), 114		
get_bath_temp()	( <i>LaudaProRp245e</i> method), 118		
get_by_p_id()	( <i>LabJack.DeviceType</i> class method), 112		
get_by_p_id()	( <i>LJMCommunicationConfig.DeviceType</i> class method), 18		
get_cal_current_source()	( <i>LabJack</i> method), 114		
get_channel_offset()	( <i>RTO1024</i> method), 152		
get_channel_position()	( <i>RTO1024</i> method), 152		
get_channel_range()	( <i>RTO1024</i> method), 152		
get_channel_scale()	( <i>RTO1024</i> method), 152		
get_channel_state()	( <i>RTO1024</i> method), 152		
get_clock()	( <i>LabJack</i> method), 114		
get_controller_information()	( <i>NewportSMC100PP</i> method), 132		
get_current()	( <i>HeinzingerDI</i> method), 105		
get_dc_volt()	( <i>ILS2T</i> method), 161		
get_device()	( <i>DeviceSequenceMixin</i> method), 73		
get_device_type()	( <i>LaudaProRp245e</i> method), 118		
get_devices()	( <i>DeviceSequenceMixin</i> method), 73		
get_digital_input()	( <i>LabJack</i> method), 114		
get_error_code()	( <i>ILS2T</i> method), 161		
get_error_queue()	( <i>VisaDevice</i> method), 173		
get_full_scale_mbar()	( <i>PfeifferTPG</i> method), 145		
get_full_scale_unitless()	( <i>PfeifferTPG</i> method), 145		
get_identification()	( <i>VisaDevice</i> method), 173		
get_interface_version()	( <i>HeinzingerDI</i> method), 105		
get_motor_configuration()	( <i>NewportSMC100PP</i> method), 132		
get_move_duration()	( <i>NewportSMC100PP</i> method), 133		

get\_negative\_software\_limit()  
*(portSMC100PP method)*, 133

get\_number\_of\_recordings()  
*(HeinzingerDI method)*, 105

get\_objects\_node()  
*(Client method)*, 21

get\_objects\_node()  
*(Server method)*, 24

get\_output()  
*(PSI9000 method)*, 86

get\_position()  
*(ILS2T method)*, 161

get\_position()  
*(NewportSMC100PP method)*, 133

get\_positive\_software\_limit()  
*(NewportSMC100PP method)*, 133

get\_product\_id()  
*(LabJack method)*, 114

get\_product\_name()  
*(LabJack method)*, 114

get\_product\_type()  
*(LabJack method)*, 115

get\_pulse\_energy\_and\_rate()  
*(CryLasLaser method)*, 79

get\_reference\_point()  
*(RTO1024 method)*, 152

get\_register()  
*(FuGProbusV method)*, 97

get\_repetitions()  
*(RTO1024 method)*, 153

get\_sbus\_rh()  
*(LabJack method)*, 115

get\_sbus\_temp()  
*(LabJack method)*, 115

get\_serial\_number()  
*(HeinzingerDI method)*, 105

get\_serial\_number()  
*(LabJack method)*, 115

get\_state()  
*(NewportSMC100PP method)*, 134

get\_status()  
*(ILS2T method)*, 161

get\_system\_lock()  
*(PSI9000 method)*, 86

get\_temperature()  
*(ILS2T method)*, 161

get\_timestamps()  
*(RTO1024 method)*, 153

get\_ui\_lower\_limits()  
*(PSI9000 method)*, 86

get\_ui\_upper\_limits()  
*(PSI9000 method)*, 86

get\_voltage()  
*(HeinzingerDI method)*, 105

get\_voltage\_current\_setpoint()  
*(PSI9000 method)*, 86

GetAttr  
*(class in hvl\_ccb.utils.conversion.utils)*, 177

go\_home()  
*(NewportSMC100PP method)*, 134

go\_to\_configuration()  
*(NewportSMC100PP method)*, 134

GREEN\_NOT\_READY  
*(SafetyStatus attribute)*, 44

GREEN\_READY  
*(SafetyStatus attribute)*, 44

## H

HARDWARE  
*(CryLasLaser.RepetitionRates attribute)*, 79

HEAD  
*(CryLasLaser.AnswersStatus attribute)*, 78

HeinzingerConfig  
*(class in hvl\_ccb.dev.heinzinger)*, 103

HeinzingerConfig.RecordingsEnum  
*(class in hvl\_ccb.dev.heinzinger)*, 103

HeinzingerDI  
*(class in hvl\_ccb.dev.heinzinger)*, 104

HeinzingerDI.OutputStatus  
*(class in hvl\_ccb.dev.heinzinger)*, 105

HeinzingerPNC  
*(class in hvl\_ccb.dev.heinzinger)*, 107

HeinzingerPNC.UnitCurrent  
*(class in hvl\_ccb.dev.heinzinger)*, 107

(New- HeinzingerPNC.UnitVoltage  
*(class in hvl\_ccb.dev.heinzinger)*, 107

HeinzingerPNCDeviceNotRecognizedError, 108

HeinzingerPNCError, 108

HeinzingerPNCMaxCurrentExceededError, 108

HeinzingerPNCMaxVoltageExceededError, 108

HeinzingerSerialCommunication  
*(class in hvl\_ccb.dev.heinzinger)*, 108

HeinzingerSerialCommunicationConfig  
*(class in hvl\_ccb.dev.heinzinger)*, 108

HIGH  
*(LabJack.DIOStatus attribute)*, 112

high\_resolution  
*(FuGProbusVSetRegisters property)*, 100

home\_search\_polling\_interval  
*(NewportSMC100PPConfig attribute)*, 138

home\_search\_timeout  
*(NewportSMC100PPConfig attribute)*, 138

home\_search\_type  
*(NewportSMC100PPConfig attribute)*, 138

home\_search\_velocity  
*(NewportSMC100PPConfig attribute)*, 138

HOME\_STARTED  
*(NewportSMC100PPSerialCommunication.ControllerError attribute)*, 140

HomeSwitch  
*(NewportSMC100PPConfig.HomeSearch attribute)*, 138

HomeSwitch\_and\_Index  
*(NewportSMC100PPConfig.HomeSearch attribute)*, 138

HOMING  
*(NewportStates attribute)*, 143

HOMING\_FROM\_RS232  
*(NewportSMC100PP.StateMessages attribute)*, 131

HOMING\_FROM\_SMC  
*(NewportSMC100PP.StateMessages attribute)*, 131

HOMING\_TIMEOUT  
*(NewportSMC100PP.MotorErrors attribute)*, 130

host  
*(ModbusTcpCommunicationConfig attribute)*, 20

host  
*(OpcUaCommunicationConfig attribute)*, 23

host  
*(TcpCommunicationConfig attribute)*, 29

host  
*(TelnetCommunicationConfig attribute)*, 31

host  
*(VisaCommunicationConfig attribute)*, 34

HT  
*(NewportConfigCommands attribute)*, 130

HUNDRED\_VOLT  
*(ACVoltageRange attribute)*, 55

HUNDRED\_VOLT  
*(DCVoltageRange attribute)*, 56

hvl\_ccb  
*module*, 185

hvl\_ccb.comm  
*module*, 35

hvl\_ccb.comm.base  
*module*, 11

hvl\_ccb.comm.labjack\_ljm  
*module*, 16

hvl\_ccb.comm.modbus\_tcp  
*module*, 19

hvl\_ccb.comm.opc  
    module, 21

hvl\_ccb.comm.serial  
    module, 24

hvl\_ccb.comm.tcp  
    module, 28

hvl\_ccb.comm.telnet  
    module, 30

hvl\_ccb.comm.visa  
    module, 32

hvl\_ccb.configuration  
    module, 181

hvl\_ccb.dev  
    module, 174

hvl\_ccb.dev.base  
    module, 72

hvl\_ccb.dev.crylas  
    module, 75

hvl\_ccb.dev.cube  
    module, 48

hvl\_ccb.dev.cube.base  
    module, 36

hvl\_ccb.dev.cube.constants  
    module, 41

hvl\_ccb.dev.cube.picube  
    module, 45

hvl\_ccb.dev.ea\_psi9000  
    module, 85

hvl\_ccb.dev.fluke884x  
    module, 57

hvl\_ccb.dev.fluke884x.base  
    module, 48

hvl\_ccb.dev.fluke884x.constants  
    module, 53

hvl\_ccb.dev.fluke884x.ranges  
    module, 54

hvl\_ccb.dev.fug  
    module, 90

hvl\_ccb.dev.heinzinger  
    module, 103

hvl\_ccb.dev.hIGHLAND\_T560  
    module, 62

hvl\_ccb.dev.hIGHLAND\_T560.base  
    module, 57

hvl\_ccb.dev.hIGHLAND\_T560.channel  
    module, 59

hvl\_ccb.dev.hIGHLAND\_T560.device  
    module, 60

hvl\_ccb.dev.labjack  
    module, 110

hvl\_ccb.dev.lauda  
    module, 118

hvl\_ccb.dev.mbw973  
    module, 125

hvl\_ccb.dev.newport  
    module, 129

hvl\_ccb.dev.pfeiffer\_tpg  
    module, 144

hvl\_ccb.dev.rs\_rto1024  
    module, 150

hvl\_ccb.dev.se\_ilS2t  
    module, 159

hvl\_ccb.dev.sst\_luminox  
    module, 166

hvl\_ccb.dev.technix  
    module, 68

hvl\_ccb.dev.technix.base  
    module, 63

hvl\_ccb.dev.technix.device  
    module, 66

hvl\_ccb.dev.utils  
    module, 171

hvl\_ccb.dev.visa  
    module, 172

hvl\_ccb.error  
    module, 183

hvl\_ccb.experiment\_manager  
    module, 183

hvl\_ccb.utils  
    module, 181

hvl\_ccb.utils.conversion  
    module, 178

hvl\_ccb.utils.conversion.map\_range  
    module, 174

hvl\_ccb.utils.conversion.sensor  
    module, 175

hvl\_ccb.utils.conversion.unit  
    module, 176

hvl\_ccb.utils.conversion.utils  
    module, 177

hvl\_ccb.utils.enum  
    module, 178

hvl\_ccb.utils.typing  
    module, 180

hvl\_ccb.utils.validation  
    module, 180

hysteresis\_compensation  
    (New-portSMC100PPConfig attribute), 138

|

ID (*FuGProbusIVCommands attribute*), 96

identification (*Fluke8845a property*), 50

Identification\_error (*PfeifferTPG.SensorStatus attribute*), 144

identifier (*LJMCommunicationConfig attribute*), 18

identify\_device() (*FuG method*), 91

identify\_device() (*HeinzingerPNC method*), 107

identify\_sensors() (*PfeifferTPG method*), 146

IKR (*PfeifferTPG.SensorTypes* attribute), 145  
 IKR11 (*PfeifferTPG.SensorTypes* attribute), 145  
 IKR9 (*PfeifferTPG.SensorTypes* attribute), 145  
 ILS2T (class in *hvl\_ccb.dev.se\_il2t*), 159  
*ILS2T.ActionsPtp* (class in *hvl\_ccb.dev.se\_il2t*), 159  
*ILS2T.Mode* (class in *hvl\_ccb.dev.se\_il2t*), 159  
*ILS2T.Ref16Jog* (class in *hvl\_ccb.dev.se\_il2t*), 159  
*ILS2T.State* (class in *hvl\_ccb.dev.se\_il2t*), 160  
*ILS2TConfig* (class in *hvl\_ccb.dev.se\_il2t*), 163  
 ILS2TError, 164  
*ILS2TModbusTcpCommunication* (class in *hvl\_ccb.dev.se\_il2t*), 164  
*ILS2TModbusTcpCommunicationConfig* (class in *hvl\_ccb.dev.se\_il2t*), 164  
*ILS2TRegAddr* (class in *hvl\_ccb.dev.se\_il2t*), 165  
*ILS2TRegDatatype* (class in *hvl\_ccb.dev.se\_il2t*), 165  
 IMMEDIATE (*TriggerSource* attribute), 54  
 IMMEDIATELY (*FuGRampModes* attribute), 101  
 IMPULSE\_140KV (*PowerSetup* attribute), 44  
 IMR (*PfeifferTPG.SensorTypes* attribute), 145  
 INACTIVE (*CryLasLaser.AnswersStatus* attribute), 78  
 INACTIVE (*DoorStatus* attribute), 42  
 INACTIVE (*EarthingStickStatus* attribute), 43  
 inhibit (*Technix* property), 66  
 init\_attenuation (*CryLasAttenuatorConfig* attribute), 76  
 init\_monitored\_nodes() (*OpcUaCommunication* method), 21  
 init\_shutter\_status (*CryLasLaserConfig* attribute), 82  
 initialize() (*NewportSMC100PP* method), 134  
 INITIALIZED (*ExperimentStatus* attribute), 185  
 INITIALIZING (*ExperimentStatus* attribute), 185  
 INITIALIZING (*SafetyStatus* attribute), 44  
 initiate\_trigger() (*Fluke8845a* method), 50  
 INPUT (*FuGProbusVRegisterGroups* attribute), 99  
 INPUT (*GateMode* attribute), 57  
 INSTALL (*AutoInstallMode* attribute), 57  
 INT32 (*ILS2TRegDatatype* attribute), 166  
 INT\_SYNTHESIZER (*TriggerMode* attribute), 59  
 interface\_type (*PSI9000VisaCommunicationConfig* attribute), 89  
 interface\_type (*RTO1024VisaCommunicationConfig* attribute), 158  
 interface\_type (*VisaCommunicationConfig* attribute), 34  
 internal (*LabJack.CjcType* attribute), 111  
 INTERNAL (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 122  
 IO\_SCANNING (*ILS2TRegAddr* attribute), 165  
 IoScanningModeValueError, 166  
 is\_configdataclass (*AsyncCommunicationProtocolConfig* attribute), 14  
 is\_configdataclass (*BaseCubeConfiguration* attribute), 39  
 is\_configdataclass (*CryLasAttenuatorConfig* attribute), 76  
 is\_configdataclass (*CryLasLaserConfig* attribute), 82  
 is\_configdataclass (*EmptyConfig* attribute), 74, 182  
 is\_configdataclass (*Fluke8845aConfig* attribute), 51  
 is\_configdataclass (*FuGConfig* attribute), 92  
 is\_configdataclass (*HeinzingerConfig* attribute), 104  
 is\_configdataclass (*ILS2TConfig* attribute), 163  
 is\_configdataclass (*LaudaProRp245eConfig* attribute), 123  
 is\_configdataclass (*LJMCommunicationConfig* attribute), 18  
 is\_configdataclass (*LuminoxConfig* attribute), 168  
 is\_configdataclass (*MBW973Config* attribute), 127  
 is\_configdataclass (*ModbusTcpCommunicationConfig* attribute), 20  
 is\_configdataclass (*NewportSMC100PPConfig* attribute), 138  
 is\_configdataclass (*OpcUaCommunicationConfig* attribute), 23  
 is\_configdataclass (*PfeifferTPGConfig* attribute), 147  
 is\_configdataclass (*T560Config* attribute), 62  
 is\_configdataclass (*TcpCommunicationConfig* attribute), 29  
 is\_configdataclass (*TechnixConfig* attribute), 68  
 is\_configdataclass (*VisaCommunicationConfig* attribute), 34  
 is\_done() (*MBW973* method), 125  
 is\_error() (*ExperimentManager* method), 184  
 is\_finished() (*ExperimentManager* method), 184  
 is\_generic\_type\_hint() (in *module hvl\_ccb.utils.typing*), 180  
 is\_in\_range() (*ILS2TRegDatatype* method), 166  
 is\_inactive (*CryLasLaser.LaserStatus* property), 79  
 is\_open (*Client* property), 21  
 is\_open (*LJMCommunication* property), 16  
 is\_open (*OpcUaCommunication* property), 22  
 is\_open (*SerialCommunication* property), 25  
 is\_open (*TelnetCommunication* property), 30  
 is\_polling() (*Poller* method), 171  
 is\_ready (*CryLasLaser.LaserStatus* property), 79  
 is\_running() (*ExperimentManager* method), 184  
 is\_started (*Technix* property), 66  
 is\_valid\_scale\_range\_reversed\_str() (*PfeifferTPGConfig.Model* method), 147

**J**

J (*LabJack.ThermocoupleType* attribute), 113  
 jerk\_time (*NewportSMC100PPConfig* attribute), 138  
 JOG (*ILS2T.Mode* attribute), 159

jog\_run() (*ILS2T method*), 162  
jog\_stop() (*ILS2T method*), 162  
JOGGING (*NewportStates attribute*), 143  
JOGGING\_FROM\_DISABLE  
    *portSMC100PP.StateMessages* (New-  
        attribute), 131  
JOGGING\_FROM\_READY  
    *portSMC100PP.StateMessages* (New-  
        attribute), 131  
JOGN\_FAST (*ILS2TRegAddr attribute*), 165  
JOGN\_SLOW (*ILS2TRegAddr attribute*), 165  
JR (*NewportConfigCommands attribute*), 130

## K

K (*LabJack.TemperatureUnit attribute*), 113  
K (*LabJack.ThermocoupleType attribute*), 113  
K (*Temperature attribute*), 177  
KELVIN (*Temperature attribute*), 177  
keys() (*AsyncCommunicationProtocolConfig class*  
    *method*), 14  
keys() (*BaseCubeConfiguration class* *method*), 39  
keys() (*BaseCubeOpcUaCommunicationConfig class*  
    *method*), 41  
keys() (*CryLasAttenuatorConfig class* *method*), 76  
keys() (*CryLasAttenuatorSerialCommunicationConfig class*  
    *method*), 77  
keys() (*CryLasLaserConfig class* *method*), 82  
keys() (*CryLasLaserSerialCommunicationConfig class*  
    *method*), 84  
keys() (*EmptyConfig class* *method*), 74, 182  
keys() (*Fluke8845aConfig class* *method*), 51  
keys() (*Fluke8845aTelnetCommunicationConfig class*  
    *method*), 52  
keys() (*FuGConfig class* *method*), 92  
keys() (*FuGSerialCommunicationConfig class* *method*),  
    102  
keys() (*HeinzingerConfig class* *method*), 104  
keys() (*HeinzingerSerialCommunicationConfig class*  
    *method*), 109  
keys() (*ILS2TConfig class* *method*), 163  
keys() (*ILS2TModbusTcpCommunicationConfig class*  
    *method*), 164  
keys() (*LaudaProRp245eConfig class* *method*), 123  
keys() (*LaudaProRp245eTcpCommunicationConfig class*  
    *method*), 124  
keys() (*LJMCommunicationConfig class* *method*), 18  
keys() (*LuminoxConfig class* *method*), 168  
keys() (*LuminoxSerialCommunicationConfig class*  
    *method*), 170  
keys() (*MBW973Config class* *method*), 127  
keys() (*MBW973SerialCommunicationConfig class*  
    *method*), 128  
keys() (*ModbusTcpCommunicationConfig class*  
    *method*), 20

**keys()** (*NewportSMC100PPConfig class method*), 138  
**keys()** (*NewportSMC100PPSerialCommunicationConfig class method*), 143  
**keys()** (*OpcUaCommunicationConfig class method*), 23  
**keys()** (*PfeifferTPGConfig class method*), 147  
**keys()** (*PfeifferTPGSerialCommunicationConfig class method*), 149  
**keys()** (*PICubeConfiguration class method*), 47  
**keys()** (*PICubeOpcUaCommunicationConfig class method*), 47  
**keys()** (*PSI9000Config class method*), 88  
**keys()** (*PSI9000VisaCommunicationConfig class method*), 90  
**keys()** (*RTO1024Config class method*), 157  
**keys()** (*RTO1024VisaCommunicationConfig class method*), 158  
**keys()** (*SerialCommunicationConfig class method*), 26  
**keys()** (*T560CommunicationConfig class method*), 58  
**keys()** (*T560Config class method*), 62  
**keys()** (*TcpCommunicationConfig class method*), 29  
**keys()** (*TechnixConfig class method*), 68  
**keys()** (*TechnixSerialCommunicationConfig class method*), 64  
**keys()** (*TechnixTelnetCommunicationConfig class method*), 65  
**keys()** (*TelnetCommunicationConfig class method*), 31  
**keys()** (*VisaCommunicationConfig class method*), 34  
**keys()** (*VisaDeviceConfig class method*), 173  
kV (*HeinzingerPNC.UnitVoltage attribute*), 107

**L**

LabJack (*class in hvl\_ccb.dev.labjack*), 111  
LabJack.AInRange (*class in hvl\_ccb.dev.labjack*), 111  
LabJack.BitLimit (*class in hvl\_ccb.dev.labjack*), 111  
LabJack.CalMicroAmpere (*class in hvl\_ccb.dev.labjack*), 111  
LabJack.CjcType (*class in hvl\_ccb.dev.labjack*), 111  
LabJack.ClockFrequency (*class in hvl\_ccb.dev.labjack*), 112  
LabJack.DeviceType (*class in hvl\_ccb.dev.labjack*), 112  
LabJack.DIOStatus (*class in hvl\_ccb.dev.labjack*), 112  
LabJack.TemperatureUnit (*class in hvl\_ccb.dev.labjack*), 113  
LabJack.ThermocoupleType (*class in hvl\_ccb.dev.labjack*), 113  
LabJackError, 117  
LabJackIdentifierDIOError, 117  
laser\_off() (*CryLasLaser method*), 80  
laser\_on() (*CryLasLaser method*), 80  
LaudaProRp245e (*class in hvl\_ccb.dev.lauda*), 118  
LaudaProRp245eCommand (*class in hvl\_ccb.dev.lauda*), 120  
LaudaProRp245eCommandError, 122

LaudaProRp245eConfig (*class in hvl\_ccb.dev.lauda*), 122  
LaudaProRp245eConfig.ExtControlModeEnum (*class in hvl\_ccb.dev.lauda*), 122  
LaudaProRp245eConfig.OperationModeEnum (*class in hvl\_ccb.dev.lauda*), 122  
LaudaProRp245eTcpCommunication (*class in hvl\_ccb.dev.lauda*), 123  
LaudaProRp245eTcpCommunicationConfig (*class in hvl\_ccb.dev.lauda*), 124  
LEM4000S (*class in hvl\_ccb.utils.conversion.sensor*), 175  
LF (*FuGTerminators attribute*), 103  
LFCR (*FuGTerminators attribute*), 103  
LINE\_1 (*MessageBoard attribute*), 43  
LINE\_10 (*MessageBoard attribute*), 43  
LINE\_11 (*MessageBoard attribute*), 43  
LINE\_12 (*MessageBoard attribute*), 43  
LINE\_13 (*MessageBoard attribute*), 43  
LINE\_14 (*MessageBoard attribute*), 43  
LINE\_15 (*MessageBoard attribute*), 43  
LINE\_2 (*MessageBoard attribute*), 43  
LINE\_3 (*MessageBoard attribute*), 43  
LINE\_4 (*MessageBoard attribute*), 43  
LINE\_5 (*MessageBoard attribute*), 43  
LINE\_6 (*MessageBoard attribute*), 43  
LINE\_7 (*MessageBoard attribute*), 43  
LINE\_8 (*MessageBoard attribute*), 43  
LINE\_9 (*MessageBoard attribute*), 43  
list\_directory() (*RTO1024 method*), 153  
LJMCommunication (*class in hvl\_ccb.comm.labjack\_ljm*), 16  
LJMCommunicationConfig (*class in hvl\_ccb.comm.labjack\_ljm*), 17  
LJMCommunicationConfig.ConnectionType (*class in hvl\_ccb.comm.labjack\_ljm*), 17  
LJMCommunicationConfig.DeviceType (*class in hvl\_ccb.comm.labjack\_ljm*), 17  
LJMCommunicationError, 18  
lm34 (*LabJack.CjcType attribute*), 112  
LMT70A (*class in hvl\_ccb.utils.conversion.sensor*), 175  
load\_configuration() (*RTO1024 method*), 153  
load\_device\_configuration() (*T560 method*), 61  
local\_display() (*RTO1024 method*), 153  
LOCKED (*DoorStatus attribute*), 42  
LOW (*LabJack.DIOStatus attribute*), 112  
LOWER\_TEMP (*LaudaProRp245eCommand attribute*), 121  
lower\_temp (*LaudaProRp245eConfig attribute*), 123  
Luminox (*class in hvl\_ccb.dev.sst\_luminox*), 167  
LuminoxConfig (*class in hvl\_ccb.dev.sst\_luminox*), 167  
LuminoxError, 168  
LuminoxMeasurementType (*class in hvl\_ccb.dev.sst\_luminox*), 168  
LuminoxMeasurementTypeDict (*in module hvl\_ccb.dev.sst\_luminox*), 169  
LuminoxMeasurementTypeError, 169  
LuminoxMeasurementTypeValue (*in module hvl\_ccb.dev.sst\_luminox*), 169  
LuminoxOutputMode (*class in hvl\_ccb.dev.sst\_luminox*), 169  
LuminoxOutputModeError, 169  
LuminoxSerialCommunication (*class in hvl\_ccb.dev.sst\_luminox*), 169  
LuminoxSerialCommunicationConfig (*class in hvl\_ccb.dev.sst\_luminox*), 170  
LUT (*LMT70A attribute*), 176

## M

mA (*HeinzingerPNC.UnitCurrent attribute*), 107  
MANUAL (*EarthingStickOperatingStatus attribute*), 42  
MapBitAsymRange (*class in hvl\_ccb.utils.conversion.map\_range*), 174  
MapBitSymRange (*class in hvl\_ccb.utils.conversion.map\_range*), 175  
MapRanges (*class in hvl\_ccb.utils.conversion.map\_range*), 175  
MARK (*SerialCommunicationParity attribute*), 27  
max\_current (*FuG property*), 91  
max\_current (*HeinzingerPNC property*), 107  
max\_current (*Technix property*), 66  
max\_current (*TechnixConfig attribute*), 68  
max\_current\_hardware (*FuG property*), 91  
max\_current\_hardware (*HeinzingerPNC property*), 107  
max\_pr\_number (*LaudaProRp245eConfig attribute*), 123  
max\_pump\_level (*LaudaProRp245eConfig attribute*), 123  
max\_timeout\_retry\_nr (*OpcUaCommunicationConfig attribute*), 23  
max\_voltage (*FuG property*), 91  
max\_voltage (*HeinzingerPNC property*), 107  
max\_voltage (*Technix property*), 67  
max\_voltage (*TechnixConfig attribute*), 68  
max\_voltage\_hardware (*FuG property*), 91  
max\_voltage\_hardware (*HeinzingerPNC property*), 107  
MAXIMUM (*LabJack.ClockFrequency attribute*), 112  
MBW973 (*class in hvl\_ccb.dev.mbw973*), 125  
MBW973Config (*class in hvl\_ccb.dev.mbw973*), 127  
MBW973ControlRunningError, 127  
MBW973Error, 127  
MBW973PumpRunningError, 128  
MBW973SerialCommunication (*class in hvl\_ccb.dev.mbw973*), 128  
MBW973SerialCommunicationConfig (*class in hvl\_ccb.dev.mbw973*), 128  
measure() (*Fluke8845a method*), 50  
measure() (*PfeifferTPG method*), 146  
measure\_all() (*PfeifferTPG method*), 146

**measure\_current()** (*HeinzingerDI method*), 106  
**measure\_voltage()** (*HeinzingerDI method*), 106  
**measure\_voltage\_current()** (*PSI9000 method*), 87  
**measurement\_function** (*Fluke8845a property*), 50  
**MeasurementFunction** (class) in *hvl\_ccb.dev.fluke884x.constants*, 53  
**MEDIUM\_FILTER** (*FilterRange attribute*), 56  
**MessageBoard** (class in *hvl\_ccb.dev(cube.constants)*), 43  
**micro\_step\_per\_full\_step\_factor** (*NewportSMC100PPConfig attribute*), 139  
**MILLIMETER\_MERCURY** (*Pressure attribute*), 176  
**MINIMUM** (*LabJack.ClockFrequency attribute*), 112  
**MMHG** (*Pressure attribute*), 176  
**ModbusTcpCommunication** (class) in *hvl\_ccb.comm.modbus\_tcp*, 19  
**ModbusTcpCommunicationConfig** (class) in *hvl\_ccb.comm.modbus\_tcp*, 20  
**ModbusTcpConnectionFailedError**, 21  
**model** (*PfeifferTPGConfig attribute*), 147  
**module**  
    *hvl\_ccb*, 185  
    *hvl\_ccb.comm*, 35  
    *hvl\_ccb.comm.base*, 11  
    *hvl\_ccb.comm.labjack\_ljm*, 16  
    *hvl\_ccb.comm.modbus\_tcp*, 19  
    *hvl\_ccb.comm.opc*, 21  
    *hvl\_ccb.comm.serial*, 24  
    *hvl\_ccb.comm.tcp*, 28  
    *hvl\_ccb.comm.telnet*, 30  
    *hvl\_ccb.comm.visa*, 32  
    *hvl\_ccb.configuration*, 181  
    *hvl\_ccb.dev*, 174  
    *hvl\_ccb.dev.base*, 72  
    *hvl\_ccb.dev.crylas*, 75  
    *hvl\_ccb.dev.cube*, 48  
    *hvl\_ccb.dev.cube.base*, 36  
    *hvl\_ccb.dev.cube.constants*, 41  
    *hvl\_ccb.dev.cube.picube*, 45  
    *hvl\_ccb.dev.ea\_psi9000*, 85  
    *hvl\_ccb.dev.fluke884x*, 57  
    *hvl\_ccb.dev.fluke884x.base*, 48  
    *hvl\_ccb.dev.fluke884x.constants*, 53  
    *hvl\_ccb.dev.fluke884x.ranges*, 54  
    *hvl\_ccb.dev.fug*, 90  
    *hvl\_ccb.dev.heinzinger*, 103  
    *hvl\_ccb.dev.highland\_t560*, 62  
    *hvl\_ccb.dev.highland\_t560.base*, 57  
    *hvl\_ccb.dev.highland\_t560.channel*, 59  
    *hvl\_ccb.dev.highland\_t560.device*, 60  
    *hvl\_ccb.dev.labjack*, 110  
    *hvl\_ccb.dev.lauda*, 118  
    *hvl\_ccb.dev.mbw973*, 125  
    *hvl\_ccb.dev.newport*, 129  
    *hvl\_ccb.dev.pfeiffer\_tpg*, 144  
  
**hvl\_ccb.dev.rs\_rto1024**, 150  
**hvl\_ccb.dev.se\_ilst2t**, 159  
**hvl\_ccb.dev.sst\_luminox**, 166  
**hvl\_ccb.dev.technix**, 68  
**hvl\_ccb.dev.technix.base**, 63  
**hvl\_ccb.dev.technix.device**, 66  
**hvl\_ccb.dev.utils**, 171  
**hvl\_ccb.dev.visa**, 172  
**hvl\_ccb.error**, 183  
**hvl\_ccb.experiment\_manager**, 183  
**hvl\_ccb.utils**, 181  
**hvl\_ccb.utils.conversion**, 178  
**hvl\_ccb.utils.conversion.map\_range**, 174  
**hvl\_ccb.utils.conversion.sensor**, 175  
**hvl\_ccb.utils.conversion.unit**, 176  
**hvl\_ccb.utils.conversion.utils**, 177  
**hvl\_ccb.utils.enum**, 178  
**hvl\_ccb.utils.typing**, 180  
**hvl\_ccb.utils.validation**, 180  
**MONITOR\_I** (*FuGProbusVRegisterGroups attribute*), 99  
**MONITOR\_V** (*FuGProbusVRegisterGroups attribute*), 99  
**most\_recent\_error** (*FuGProbusVConfigRegisters property*), 97  
**motion\_distance\_per\_full\_step** (*NewportSMC100PPConfig attribute*), 139  
**motor\_config** (*NewportSMC100PPConfig property*), 139  
**move\_to\_absolute\_position()** (*NewportSMC100PP method*), 135  
**move\_to\_relative\_position()** (*NewportSMC100PP method*), 135  
**move\_wait\_sec** (*NewportSMC100PPConfig attribute*), 139  
**MOVING** (*NewportSMC100PP.StateMessages attribute*), 131  
**MOVING** (*NewportStates attribute*), 143  
**MS\_NOMINAL\_CURRENT** (*PSI9000 attribute*), 86  
**MS\_NOMINAL\_VOLTAGE** (*PSI9000 attribute*), 86  
**MULTI\_COMMANDS\_MAX** (*VisaCommunication attribute*), 32  
**MULTI\_COMMANDS\_SEPARATOR** (*VisaCommunication attribute*), 32

## N

**name** (*Fluke8845aConfig attribute*), 51  
**NameEnum** (class in *hvl\_ccb.utils.enum*), 178  
**NAMES** (*SerialCommunicationParity attribute*), 27  
**names()** (*RTO1024.TriggerModes class method*), 151  
**namespace\_index** (*BaseCubeConfiguration attribute*), 39  
**NED\_END\_OF\_TURN** (*NewportSMC100PP.MotorErrors attribute*), 130  
**NEG** (*ILS2T.Ref16Jog attribute*), 160  
**NEG\_FAST** (*ILS2T.Ref16Jog attribute*), 160

NEGATIVE ( <i>FuGPolarities attribute</i> ), 95			
NEGATIVE ( <i>Polarity attribute</i> ), 44			
negative_software_limit	(New-	NO_REF_FROM_RESET	(New-
<i>portSMC100PPConfig attribute</i> ), 139	portSMC100PP.StateMessages	portSMC100PP.StateMessages	attribute),
NewportConfigCommands	(class	in	131
<i>hvl_ccb.dev.newport</i> ), 129			
NewportControllerError	, 130	No_sensor ( <i>PfeifferTPG.SensorStatus attribute</i> ), 144	
NewportError	, 130	NO_SOURCE ( <i>PowerSetup attribute</i> ), 44	
NewportMotorError	, 130	noise_level_measurement_channel_1 ( <i>BaseCube-</i>	
NewportMotorPowerSupplyWasCutError	, 130	<i>Configuration attribute</i> ), 40	
NewportSerialCommunicationError	, 143	noise_level_measurement_channel_2 ( <i>BaseCube-</i>	
NewportSMC100PP ( <i>class in hvl_ccb.dev.newport</i> ), 130		<i>Configuration attribute</i> ), 40	
NewportSMC100PP.MotorErrors	(class	noise_level_measurement_channel_3 ( <i>BaseCube-</i>	
<i>hvl_ccb.dev.newport</i> ), 130	in	<i>Configuration attribute</i> ), 40	
NewportSMC100PP.StateMessages	(class	noise_level_measurement_channel_4 ( <i>BaseCube-</i>	
<i>hvl_ccb.dev.newport</i> ), 131	in	<i>Configuration attribute</i> ), 40	
NewportSMC100PPConfig	(class	NONE ( <i>ILS2T.Ref16Jog attribute</i> ), 160	
<i>hvl_ccb.dev.newport</i> ), 137	in	NONE ( <i>LabJack.ThermocoupleType attribute</i> ), 113	
NewportSMC100PPConfig.EspStageConfig	(class	None ( <i>PfeifferTPG.SensorTypes attribute</i> ), 145	
<i>hvl_ccb.dev.newport</i> ), 137	in	NONE ( <i>SerialCommunicationParity attribute</i> ), 27	
NewportSMC100PPConfig.HomeSearch	(class	NORMAL ( <i>RTO1024.TriggerModes attribute</i> ), 151	
<i>hvl_ccb.dev.newport</i> ), 138	in	noSen ( <i>PfeifferTPG.SensorTypes attribute</i> ), 145	
NewportSMC100PPSerialCommunication	(class	noSENSOR ( <i>PfeifferTPG.SensorTypes attribute</i> ), 145	
<i>hvl_ccb.dev.newport</i> ), 139	in	nr_trials_activate ( <i>LuminoxConfig attribute</i> ), 168	
NewportSMC100PPSerialCommunication.Controller		NullCommunicationProtocol	(class
<i>(class in hvl_ccb.dev.newport)</i> , 139		<i>hvl_ccb.comm.base</i> ), 14	in
NewportSMC100PPSerialCommunicationConfig		Number	(in module <i>hvl_ccb.utils.typing</i> ), 180
<i>(class in hvl_ccb.dev.newport)</i> , 142		number_of_decimals ( <i>HeinzingerConfig attribute</i> ), 104	
NewportStates	(class	number_of_sensors ( <i>PfeifferTPG property</i> ), 146	
<i>in hvl_ccb.dev.newport</i> ), 143			
NewportUncertainPositionError	, 144		
NO ( <i>FuDigitalVal attribute</i> ), 93			
NO_ERROR ( <i>NewportSMC100PPSerialCommunication.Controller</i>			
<i>attribute</i> ), 140			
NO_REF ( <i>NewportStates attribute</i> ), 143		O	
NO_REF_ESP_STAGE_ERROR		ODD ( <i>SerialCommunicationParity attribute</i> ), 27	
<i>portSMC100PP.StateMessages</i>		OFF ( <i>AutoInstallMode attribute</i> ), 57	
131	(New-	OFF ( <i>FuDigitalVal attribute</i> ), 93	
NO_REF_FROM_CONFIG		OFF ( <i>GateMode attribute</i> ), 57	
<i>portSMC100PP.StateMessages</i>		OFF ( <i>HeinzingerDI.OutputStatus attribute</i> ), 105	
131	(New-	OFF ( <i>TriggerMode attribute</i> ), 59	
NO_REF_FROM_DISABLED		OH ( <i>NewportConfigCommands attribute</i> ), 130	
<i>portSMC100PP.StateMessages</i>		Ok ( <i>PfeifferTPG.SensorStatus attribute</i> ), 144	
131	(New-	on ( <i>FuG property</i> ), 91	
NO_REF_FROM_HOMING		ON ( <i>FuDigitalVal attribute</i> ), 93	
<i>portSMC100PP.StateMessages</i>		on ( <i>FuGProbusVDIRegisters property</i> ), 98	
131	(New-	ON ( <i>HeinzingerDI.OutputStatus attribute</i> ), 105	
NO_REF_FROM_JOGGING		ON ( <i>ILS2T.State attribute</i> ), 160	
<i>portSMC100PP.StateMessages</i>		ONE ( <i>HeinzingerConfig.RecordingsEnum attribute</i> ), 104	
131	(New-	ONE ( <i>LabJack.AInRange attribute</i> ), 111	
NO_REF_FROM_MOVING		ONE ( <i>SerialCommunicationStopbits attribute</i> ), 27	
<i>portSMC100PP.StateMessages</i>		ONE_AMPERE ( <i>ACCurrentRange attribute</i> ), 54	
131	(New-	ONE_AMPERE ( <i>DCCurrentRange attribute</i> ), 55	
NO_REF_FROM_READY		ONE_HUNDRED_MICRO_AMPERE	( <i>DCCurrentRange</i>
<i>portSMC100PP.StateMessages</i>		attribute), 55	
	(New-	ONE_HUNDRED_MILLI_AMPERE	( <i>ACCurrentRange</i>
	attribute),	attribute), 54	
		ONE_HUNDRED_MILLI_AMPERE	( <i>DCCurrentRange</i>
		attribute), 55	

ONE\_HUNDRED\_MILLI\_SECOND (*ApertureRange attribute*), 55  
ONE\_HUNDRED\_MILLI\_VOLT (*ACVoltageRange attribute*), 55  
ONE\_HUNDRED\_MILLI\_VOLT (*DCVoltageRange attribute*), 56  
ONE\_HUNDRED\_MILLION\_OHM (*ResistanceRange attribute*), 56  
ONE\_HUNDRED\_OHM (*ResistanceRange attribute*), 56  
ONE\_HUNDRED\_THOUSAND\_OHM (*ResistanceRange attribute*), 56  
ONE\_HUNDREDTH (*LabJack.AInRange attribute*), 111  
ONE\_MILLI\_AMPERE (*DCCurrentRange attribute*), 55  
ONE\_MILLION\_OHM (*ResistanceRange attribute*), 56  
ONE\_POINT\_FIVE (*SerialCommunicationStopbits attribute*), 27  
ONE\_SECOND (*ApertureRange attribute*), 55  
ONE\_TENTH (*LabJack.AInRange attribute*), 111  
ONE\_THOUSAND\_OHM (*ResistanceRange attribute*), 56  
ONE\_THOUSAND\_VOLT (*DCVoltageRange attribute*), 56  
ONE\_VOLT (*ACVoltageRange attribute*), 55  
ONE\_VOLT (*DCVoltageRange attribute*), 56  
ONLYUPWARDSOFFTOZERO (*FuGRampModes attribute*), 101  
OPC\_MAX\_YEAR (*BaseCube attribute*), 36  
OPC\_MIN\_YEAR (*BaseCube attribute*), 36  
OpcUaCommunication (*class in hvl\_ccb.comm.opc*), 21  
OpcUaCommunicationConfig (*class in hvl\_ccb.comm.opc*), 22  
OpcUaCommunicationIOError, 24  
OpcUaCommunicationTimeoutError, 24  
OpcUaSubHandler (*class in hvl\_ccb.comm.opc*), 24  
OPEN (*DoorStatus attribute*), 42  
OPEN (*EarthingStickOperation attribute*), 43  
OPEN (*EarthingStickStatus attribute*), 43  
open() (*CommunicationProtocol method*), 14  
open() (*LaudaProRp245eTcpCommunication method*), 124  
open() (*LJMCommunication method*), 16  
open() (*ModbusTcpCommunication method*), 19  
open() (*NullCommunicationProtocol method*), 15  
open() (*OpcUaCommunication method*), 22  
open() (*SerialCommunication method*), 25  
open() (*Tcp method*), 28  
open() (*TelnetCommunication method*), 30  
open() (*VisaCommunication method*), 33  
open\_interlock (*Technix property*), 67  
open\_shutter() (*CryLasLaser method*), 80  
open\_timeout (*VisaCommunicationConfig attribute*), 34  
OPENED (*CryLasLaser.AnswersShutter attribute*), 78  
OPENED (*CryLasLaserShutterStatus attribute*), 85  
operate (*BaseCube property*), 37  
operate (*PICube property*), 46  
OPERATION\_MODE (*LaudaProRp245eCommand attribute*), 121  
operation\_mode (*LaudaProRp245eConfig attribute*), 123  
optional\_defaults() (*AsyncCommunicationProtocolConfig class method*), 14  
optional\_defaults() (*BaseCubeConfiguration class method*), 40  
optional\_defaults() (*BaseCubeOpcUaCommunicationConfig class method*), 41  
optional\_defaults() (*CryLasAttenuatorConfig class method*), 76  
optional\_defaults() (*CryLasAttenuatorSerialCommunicationConfig class method*), 78  
optional\_defaults() (*CryLasLaserConfig class method*), 82  
optional\_defaults() (*CryLasLaserSerialCommunicationConfig class method*), 84  
optional\_defaults() (*EmptyConfig class method*), 74, 182  
optional\_defaults() (*Fluke8845aConfig class method*), 51  
optional\_defaults() (*Fluke8845aTelnetCommunicationConfig class method*), 52  
optional\_defaults() (*FuGConfig class method*), 93  
optional\_defaults() (*FuGSerialCommunicationConfig class method*), 102  
optional\_defaults() (*HeinzingerConfig class method*), 104  
optional\_defaults() (*HeinzingerSerialCommunicationConfig class method*), 109  
optional\_defaults() (*ILS2TConfig class method*), 163  
optional\_defaults() (*ILS2TModbusTcpCommunicationConfig class method*), 165  
optional\_defaults() (*LaudaProRp245eConfig class method*), 123  
optional\_defaults() (*LaudaProRp245eTcpCommunicationConfig class method*), 124  
optional\_defaults() (*LJMCommunicationConfig class method*), 18  
optional\_defaults() (*LuminoxConfig class method*), 168  
optional\_defaults() (*LuminoxSerialCommunicationConfig class method*), 170  
optional\_defaults() (*MBW973Config class method*), 127  
optional\_defaults() (*MBW973SerialCommunicationConfig class method*), 128  
optional\_defaults() (*ModbusTcpCommunication-*

*Config class method), 20*  
**optional\_defaults()** (*NewportSMC100PPConfig class method*), 139  
**optional\_defaults()** (*NewportSMC100PPSerialCommunicationConfig class method*), 143  
**optional\_defaults()** (*OpcUaCommunicationConfig class method*), 23  
**optional\_defaults()** (*PfeifferTPGConfig class method*), 148  
**optional\_defaults()** (*PfeifferTPGSerialCommunicationConfig class method*), 149  
**optional\_defaults()** (*PICubeConfiguration class method*), 47  
**optional\_defaults()** (*PICubeOpcUaCommunicationConfig class method*), 48  
**optional\_defaults()** (*PSI9000Config class method*), 88  
**optional\_defaults()** (*PSI9000VisaCommunicationConfig class method*), 90  
**optional\_defaults()** (*RTO1024Config class method*), 157  
**optional\_defaults()** (*RTO1024VisaCommunicationConfig class method*), 158  
**optional\_defaults()** (*SerialCommunicationConfig class method*), 26  
**optional\_defaults()** (*T560CommunicationConfig class method*), 58  
**optional\_defaults()** (*T560Config class method*), 62  
**optional\_defaults()** (*TcpCommunicationConfig class method*), 29  
**optional\_defaults()** (*TechnixConfig class method*), 68  
**optional\_defaults()** (*TechnixSerialCommunicationConfig class method*), 64  
**optional\_defaults()** (*TechnixTelnetCommunicationConfig class method*), 65  
**optional\_defaults()** (*TelnetCommunicationConfig class method*), 31  
**optional\_defaults()** (*VisaCommunicationConfig class method*), 34  
**optional\_defaults()** (*VisaDeviceConfig class method*), 174  
**OT** (*NewportConfigCommands attribute*), 130  
**out** (*FuGProbusVDORegisters property*), 99  
**OUTPUT** (*FuGProbusIVCommands attribute*), 96  
**OUTPUT** (*GateMode attribute*), 57  
**output** (*Technix property*), 67  
**output\_off()** (*FuGProbusIV method*), 96  
**output\_off()** (*HeinzingerDI method*), 106  
**output\_on()** (*HeinzingerDI method*), 106  
**OUTPUT\_POWER\_EXCEEDED** (*NewportSMC100PP.MotorErrors attribute*), 131  
**output\_status** (*HeinzingerDI property*), 106  
**OUTPUTONCMD** (*FuGProbusVRegisterGroups attribute*), 99  
**OUTPUTX0** (*FuGProbusVRegisterGroups attribute*), 99  
**OUTPUTX1** (*FuGProbusVRegisterGroups attribute*), 100  
**OUTPUTX2** (*FuGProbusVRegisterGroups attribute*), 100  
**OUTPUTXCMD** (*FuGProbusVRegisterGroups attribute*), 100  
**outX0** (*FuG property*), 91  
**outX1** (*FuG property*), 91  
**outX2** (*FuG property*), 92  
**outXCMD** (*FuG property*), 92  
**Overrange** (*PfeifferTPG.SensorStatus attribute*), 144

## P

**PA** (*Pressure attribute*), 176  
**PARAM\_MISSING\_OR\_INVALID** (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 140  
**parity** (*CryLasAttenuatorSerialCommunicationConfig attribute*), 78  
**parity** (*CryLasLaserSerialCommunicationConfig attribute*), 85  
**parity** (*FuGSerialCommunicationConfig attribute*), 102  
**parity** (*HeinzigerSerialCommunicationConfig attribute*), 109  
**parity** (*LuminoxSerialCommunicationConfig attribute*), 170  
**parity** (*MBW973SerialCommunicationConfig attribute*), 129  
**parity** (*NewportSMC100PPSerialCommunicationConfig attribute*), 143  
**parity** (*PfeifferTPGSerialCommunicationConfig attribute*), 149  
**Parity** (*SerialCommunicationConfig attribute*), 26  
**parity** (*SerialCommunicationConfig attribute*), 27  
**parse\_read\_measurement\_value()** (*LuminoxMeasurementType method*), 169  
**partial\_pressure\_o2** (*LuminoxMeasurementType attribute*), 169  
**PASCAL** (*Pressure attribute*), 176  
**pause()** (*LaudaProRp245e method*), 118  
**pause\_ramp()** (*LaudaProRp245e method*), 118  
**PBR** (*PfeifferTPG.SensorTypes attribute*), 145  
**PEAK\_CURRENT\_LIMIT** (*NewportSMC100PP.MotorErrors attribute*), 131  
**peak\_output\_current\_limit** (*NewportSMC100PPConfig attribute*), 139  
**percent\_o2** (*LuminoxMeasurementType attribute*), 169  
**PERIOD** (*MeasurementFunction attribute*), 53  
**period** (*T560 property*), 61

period_aperture ( <i>Fluke8845a attribute</i> ), 50				
PfeifferTPG ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 144				
PfeifferTPG.SensorStatus ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 144	<i>in</i>		POSITIVE ( <i>FuGPolarities attribute</i> ), 95	
PfeifferTPG.SensorTypes ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 145	<i>in</i>		POSITIVE ( <i>Polarity attribute</i> ), 44	
PfeifferTPGConfig ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 147			positive_software_limit ( <i>NewportSMC100PPConfig attribute</i> ), 139	(NewportSMC100PPConfig method), 139
PfeifferTPGConfig.Model ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 147	<i>in</i>		post_stop_pause_sec ( <i>TechnixConfig attribute</i> ), 68	
PfeifferTPGError, 148			POUNDS_PER_SQUARE_INCH ( <i>Pressure attribute</i> ), 176	
PfeifferTPGSerialCommunication ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 148	<i>in</i>		POWER_INVERTER_220V ( <i>PowerSetup attribute</i> ), 44	
PfeifferTPGSerialCommunicationConfig ( <i>class in hvl_ccb.dev.pfeiffer_tpg</i> ), 148			power_limit ( <i>PSI9000Config attribute</i> ), 88	
PICube ( <i>class in hvl_ccb.dev.cube.picube</i> ), 45			power_setup ( <i>PICube property</i> ), 46	
PICubeConfiguration ( <i>class in hvl_ccb.dev.cube.picube</i> ), 46	<i>in</i>		PowerSetup ( <i>class in hvl_ccb.dev.cube.constants</i> ), 44	
PICubeOpcUaCommunication ( <i>class in hvl_ccb.dev.cube.picube</i> ), 47	<i>in</i>		prepare_ultra_segmentation() ( <i>RTO1024 method</i> ), 153	
PICubeOpcUaCommunicationConfig ( <i>class in hvl_ccb.dev.cube.picube</i> ), 47	<i>in</i>		preserve_type() ( <i>in module hvl_ccb.utils.conversion.utils</i> ), 177	
PICubeTestParameterError, 43			Pressure ( <i>class in hvl_ccb.utils.conversion.unit</i> ), 176	
PKR ( <i>PfeifferTPG.SensorTypes attribute</i> ), 145			PSI ( <i>Pressure attribute</i> ), 177	
Polarity ( <i>class in hvl_ccb.dev.cube.constants</i> ), 43			PSI9000 ( <i>class in hvl_ccb.dev.ea_psi9000</i> ), 85	
Polarity ( <i>class in hvl_ccb.dev.hIGHLAND_T560.base</i> ), 57			PSI9000Config ( <i>class in hvl_ccb.dev.ea_psi9000</i> ), 88	
POLARITY ( <i>FuGProbusIVCommands attribute</i> ), 96			PSI9000Error, 89	
polarity ( <i>PICube property</i> ), 46			PSI9000VisaCommunication ( <i>class in hvl_ccb.dev.ea_psi9000</i> ), 89	
Poller ( <i>class in hvl_ccb.dev.utils</i> ), 171			PSI9000VisaCommunicationConfig ( <i>class in hvl_ccb.dev.ea_psi9000</i> ), 89	
polling ( <i>LuminoxOutputMode attribute</i> ), 169			PT100 ( <i>LabJack.ThermocoupleType attribute</i> ), 113	
polling_delay_sec ( <i>BaseCubeConfiguration attribute</i> ), 40	<i>at-</i>		PT1000 ( <i>LabJack.ThermocoupleType attribute</i> ), 113	
polling_interval ( <i>MBW973Config attribute</i> ), 127			PT500 ( <i>LabJack.ThermocoupleType attribute</i> ), 113	
polling_interval_sec ( <i>BaseCubeConfiguration attribute</i> ), 40			PTP ( <i>ILS2T.Mode attribute</i> ), 159	
polling_interval_sec ( <i>TechnixConfig attribute</i> ), 68			pump_init ( <i>LaudaProRp245eConfig attribute</i> ), 123	
polling_period ( <i>CryLasLaserConfig attribute</i> ), 82			PUMP_LEVEL ( <i>LaudaProRp245eCommand attribute</i> ), 121	
polling_timeout ( <i>CryLasLaserConfig attribute</i> ), 82				
port ( <i>Fluke8845aTelnetCommunicationConfig attribute</i> ), 52				
port ( <i>ModbusTcpCommunicationConfig attribute</i> ), 20				
port ( <i>OpcUaCommunicationConfig attribute</i> ), 23				
port ( <i>SerialCommunicationConfig attribute</i> ), 27				
port ( <i>T560CommunicationConfig attribute</i> ), 59				
port ( <i>TcpCommunicationConfig attribute</i> ), 29				
port ( <i>TechnixTelnetCommunicationConfig attribute</i> ), 65				
port ( <i>TelnetCommunicationConfig attribute</i> ), 31				
port ( <i>VisaCommunicationConfig attribute</i> ), 34				
POS ( <i>ILS2T.Ref16Jog attribute</i> ), 160				
POS_END_OF_TURN ( <i>NewportSMC100PP.MotorErrors attribute</i> ), 131				
POS_FAST ( <i>ILS2T.Ref16Jog attribute</i> ), 160				
POSITION ( <i>ILS2TRegAddr attribute</i> ), 165				
POSITION_OUT_OF_LIMIT ( <i>NewportSMC100PPSerialCommunication.ControllerErrors attribute</i> ), 140				

## Q

QIL ( <i>NewportConfigCommands attribute</i> ), 130				
QUERY ( <i>FuGProbusIVCommands attribute</i> ), 96				
query() ( <i>CryLasLaserSerialCommunication method</i> ), 83				
query() ( <i>Fluke8845aTelnetCommunication method</i> ), 52				
query() ( <i>FuGSerialCommunication method</i> ), 101				
query() ( <i>NewportSMC100PPSerialCommunication method</i> ), 140				
query() ( <i>PfeifferTPGSerialCommunication method</i> ), 148				
query() ( <i>SyncCommunicationProtocol method</i> ), 15				
query() ( <i>T560Communication method</i> ), 58				
query() ( <i>VisaCommunication method</i> ), 33				
query_all() ( <i>CryLasLaserSerialCommunication method</i> ), 83				
query_command() ( <i>LaudaProRp245eTcpCommunication method</i> ), 124				

**query\_multiple()** (*NewportSMC100PPSerialCommunication method*), 141  
**query\_polling()** (*Luminox method*), 167  
**query\_status()** (*Technix method*), 67  
**QUEUE** (*AutoInstallMode attribute*), 57  
**QUICK\_STOP** (*SafetyStatus attribute*), 44  
**QUICKSTOP** (*ILS2T.State attribute*), 160  
**quickstop()** (*ILS2T method*), 162  
**quit\_error()** (*BaseCube method*), 38

**R**

**R** (*LabJack.ThermocoupleType attribute*), 113  
**raise\_()** (*FuGErrorcodes method*), 95  
**RAMP\_ACC** (*ILS2TRegAddr attribute*), 165  
**RAMP\_CONTINUE** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_DECEL** (*ILS2TRegAddr attribute*), 165  
**RAMP\_DELETE** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_ITERATIONS** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_N\_MAX** (*ILS2TRegAddr attribute*), 165  
**RAMP\_PAUSE** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_SELECT** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_SET** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_START** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_STOP** (*LaudaProRp245eCommand attribute*), 121  
**RAMP\_TYPE** (*ILS2TRegAddr attribute*), 165  
**rampmode** (*FuGProbusVSetRegisters property*), 100  
**ramprate** (*FuGProbusVSetRegisters property*), 100  
**rampstate** (*FuGProbusVSetRegisters property*), 100  
**RAMPUPWARDS** (*FuGRampModes attribute*), 101  
**RangeEnum** (*class in hvl\_ccb.utils.enum*), 179  
**RATEDCURRENT** (*FuGReadbackChannels attribute*), 101  
**RATEDVOLTAGE** (*FuGReadbackChannels attribute*), 101  
**read()** (*AsyncCommunicationProtocol method*), 11  
**read()** (*BaseCube method*), 38  
**read()** (*CryLasLaserSerialCommunication method*), 84  
**read()** (*LaudaProRp245eTcpCommunication method*), 124  
**read()** (*MBW973 method*), 126  
**read()** (*OpcUaCommunication method*), 22  
**read()** (*Tcp method*), 28  
**read\_all()** (*AsyncCommunicationProtocol method*), 12  
**read\_bytes()** (*AsyncCommunicationProtocol method*), 12  
**read\_bytes()** (*SerialCommunication method*), 25  
**read\_bytes()** (*TelnetCommunication method*), 30  
**read\_float()** (*MBW973 method*), 126  
**read\_holding\_registers()** (*ModbusTcpCommunication method*), 19  
**read\_input\_registers()** (*ModbusTcpCommunication method*), 19  
**read\_int()** (*MBW973 method*), 126  
**read\_measurement()** (*RTO1024 method*), 154  
**read\_measurements()** (*MBW973 method*), 126  
**read\_name()** (*LJMCommunication method*), 16  
**read\_nonempty()** (*AsyncCommunicationProtocol method*), 12  
**read\_output\_while\_polling** (*TechnixConfig attribute*), 68  
**read\_resistance()** (*LabJack method*), 115  
**read\_single\_bytes()** (*SerialCommunication method*), 25  
**read\_streaming()** (*Luminox method*), 167  
**read\_termination** (*VisaCommunicationConfig attribute*), 35  
**read\_text()** (*AsyncCommunicationProtocol method*), 12  
**read\_text()** (*NewportSMC100PPSerialCommunication method*), 141  
**read\_text\_nonempty()** (*AsyncCommunicationProtocol method*), 12  
**READ\_TEXT\_SKIP\_PREFIXES** (*CryLasLaserSerialCommunication attribute*), 83  
**read\_thermocouple()** (*LabJack method*), 115  
**readback\_data** (*FuGProbusVConfigRegisters property*), 97  
**READBACKCHANNEL** (*FuGProbusIVCommands attribute*), 96  
**ready** (*BaseCube property*), 38  
**READY** (*CryLasLaser.AnswersStatus attribute*), 78  
**READY** (*ILS2T.State attribute*), 160  
**READY** (*NewportStates attribute*), 144  
**READY\_ACTIVE** (*CryLasLaser.LaserStatus attribute*), 79  
**READY\_FROM\_DISABLE** (*NewportSMC100PP.StateMessages attribute*), 131  
**READY\_FROM\_HOMING** (*NewportSMC100PP.StateMessages attribute*), 131  
**READY\_FROM\_JOGGING** (*NewportSMC100PP.StateMessages attribute*), 131  
**READY\_FROM\_MOVING** (*NewportSMC100PP.StateMessages attribute*), 131  
**READY\_INACTIVE** (*CryLasLaser.LaserStatus attribute*), 79  
**RED\_OPERATE** (*SafetyStatus attribute*), 44  
**RED\_READY** (*SafetyStatus attribute*), 45  
**reg\_3** (*FuGProbusVDIRegisters property*), 98  
**RegAddr** (*ILS2T attribute*), 160  
**RegDatatype** (*ILS2T attribute*), 160  
**register\_pulse\_time** (*TechnixConfig attribute*), 68

RELATIVE\_POSITION\_MOTOR (*ILS2T.ActionsPtp* attribute), 159  
RELATIVE\_POSITION\_TARGET (*ILS2T.ActionsPtp* attribute), 159  
remote (*Technix* property), 67  
remove\_device() (*DeviceSequenceMixin* method), 73  
required\_keys() (*AsyncCommunicationProtocolConfig* class method), 14  
required\_keys() (*BaseCubeConfiguration* class method), 40  
required\_keys() (*BaseCubeOpcUaCommunicationConfig* class method), 41  
required\_keys() (*CryLasAttenuatorConfig* class method), 76  
required\_keys() (*CryLasAttenuatorSerialCommunicationConfig* class method), 78  
required\_keys() (*CryLasLaserConfig* class method), 82  
required\_keys() (*CryLasLaserSerialCommunicationConfig* class method), 85  
required\_keys() (*EmptyConfig* class method), 74, 182  
required\_keys() (*Fluke8845aConfig* class method), 51  
required\_keys() (*Fluke8845aTelnetCommunicationConfig* class method), 52  
required\_keys() (*FuGConfig* class method), 93  
required\_keys() (*FuGSerialCommunicationConfig* class method), 102  
required\_keys() (*HeinzingerConfig* class method), 104  
required\_keys() (*HeinzingerSerialCommunicationConfig* class method), 109  
required\_keys() (*ILS2TConfig* class method), 164  
required\_keys() (*ILS2TModbusTcpCommunicationConfig* class method), 165  
required\_keys() (*LaudaProRp245eConfig* class method), 123  
required\_keys() (*LaudaProRp245eTcpCommunicationConfig* class method), 124  
required\_keys() (*LJMCommunicationConfig* class method), 18  
required\_keys() (*LuminoxConfig* class method), 168  
required\_keys() (*LuminoxSerialCommunicationConfig* class method), 170  
required\_keys() (*MBW973Config* class method), 127  
required\_keys() (*MBW973SerialCommunicationConfig* class method), 129  
required\_keys() (*ModbusTcpCommunicationConfig* class method), 20  
required\_keys() (*NewportSMC100PPConfig* class method), 139  
required\_keys() (*NewportSMC100PPSerialCommunicationConfig* class method), 143  
required\_keys() (*OpcUaCommunicationConfig* class method), 23  
required\_keys() (*PfeifferTPGConfig* class method), 148  
required\_keys() (*PfeifferTPGSerialCommunicationConfig* class method), 149  
required\_keys() (*PICubeConfiguration* class method), 47  
required\_keys() (*PICubeOpcUaCommunicationConfig* class method), 48  
required\_keys() (*PSI9000Config* class method), 89  
required\_keys() (*PSI9000VisaCommunicationConfig* class method), 90  
required\_keys() (*RTO1024Config* class method), 158  
required\_keys() (*RTO1024VisaCommunicationConfig* class method), 158  
required\_keys() (*SerialCommunicationConfig* class method), 27  
required\_keys() (*T560CommunicationConfig* class method), 59  
required\_keys() (*T560Config* class method), 62  
required\_keys() (*TcpCommunicationConfig* class method), 29  
required\_keys() (*TechnixConfig* class method), 68  
required\_keys() (*TechnixSerialCommunicationConfig* class method), 64  
required\_keys() (*TechnixTelnetCommunicationConfig* class method), 65  
required\_keys() (*TelnetCommunicationConfig* class method), 31  
required\_keys() (*VisaCommunicationConfig* class method), 35  
required\_keys() (*VisaDeviceConfig* class method), 174  
RESET (*FuGProbusIVCommands* attribute), 96  
reset() (*Fluke8845a* method), 50  
reset() (*FuGProbusIV* method), 96  
reset() (*NewportSMC100PP* method), 135  
reset() (*VisaDevice* method), 173  
reset\_error() (*ILS2T* method), 162  
reset\_interface() (*HeinzingerDI* method), 106  
reset\_ramp() (*LaudaProRp245e* method), 118  
ResistanceRange (class in *hvl\_ccb.dev.fluke884x.ranges*), 56  
response\_sleep\_time (*CryLasAttenuatorConfig* attribute), 77  
RMS\_CURRENT\_LIMIT (*NewportSMC100PP.MotorErrors* attribute), 131  
rpm\_max\_init (*ILS2TConfig* attribute), 164  
rs485\_address (*NewportSMC100PPConfig* attribute), 139  
RT01024 (class in *hvl\_ccb.dev.rs\_rto1024*), 150  
RT01024.TriggerModes (class in *hvl\_ccb.dev.rs\_rto1024*), 151

RT01024Config ( <i>class in hvl_ccb.dev.rs_rto1024</i> ), 157		set_acceleration() ( <i>NewportSMC100PP method</i> ), 135
RT01024Error, 158		set_acquire_length() ( <i>RTO1024 method</i> ), 154
RT01024VisaCommunication ( <i>class in hvl_ccb.dev.rs_rto1024</i> ), 158		set_ain_differential() ( <i>LabJack method</i> ), 116
RT01024VisaCommunicationConfig ( <i>class in hvl_ccb.dev.rs_rto1024</i> ), 158		set_ain_range() ( <i>LabJack method</i> ), 116
run() ( <i>ExperimentManager method</i> ), 184		set_ain_resistance() ( <i>LabJack method</i> ), 116
run() ( <i>LaudaProRp245e method</i> ), 119		set_ain_resolution() ( <i>LabJack method</i> ), 116
run_continuous_acquisition() ( <i>RTO1024 method</i> ), 154		set_ain_thermocouple() ( <i>LabJack method</i> ), 116
run_single_acquisition() ( <i>RTO1024 method</i> ), 154		set_analog_output() ( <i>LabJack method</i> ), 117
RUNNING ( <i>ExperimentStatus attribute</i> ), 185		set_attenuation() ( <i>CryLasAttenuator method</i> ), 75
<b>S</b>		set_channel_offset() ( <i>RTO1024 method</i> ), 154
S ( <i>LabJack.ThermocoupleType attribute</i> ), 113		set_channel_position() ( <i>RTO1024 method</i> ), 155
SA ( <i>NewportConfigCommands attribute</i> ), 130		set_channel_range() ( <i>RTO1024 method</i> ), 155
SafetyStatus ( <i>class in hvl_ccb.dev.cube.constants</i> ), 44		set_channel_scale() ( <i>RTO1024 method</i> ), 155
save_configuration() ( <i>RTO1024 method</i> ), 154		set_channel_state() ( <i>RTO1024 method</i> ), 156
save_device_configuration() ( <i>T560 method</i> ), 61		set_clock() ( <i>LabJack method</i> ), 117
save_waveform_history() ( <i>RTO1024 method</i> ), 154		set_control_mode() ( <i>LaudaProRp245e method</i> ), 119
SCALE ( <i>ILS2TRegAddr attribute</i> ), 165		set_current() ( <i>HeinzingerDI method</i> ), 106
ScalingFactorValueError, 166		set_current() ( <i>HeinzingerPNC method</i> ), 108
screw_scaling ( <i>NewportSMC100PPConfig attribute</i> ), 139		set_digital_output() ( <i>LabJack method</i> ), 117
send_command() ( <i>NewportSMC100PPSerialCommunication method</i> ), 141		set_external_temp() ( <i>LaudaProRp245e method</i> ), 119
send_command() ( <i>PfeifferTPGSerialCommunication method</i> ), 148		set_full_scale_mbar() ( <i>PfeifferTPG method</i> ), 146
send_hello() ( <i>Client method</i> ), 21		set_full_scale_unitless() ( <i>PfeifferTPG method</i> ), 146
send_pulses() ( <i>LabJack method</i> ), 115		set_init_attenuation() ( <i>CryLasAttenuator method</i> ), 75
send_stop() ( <i>NewportSMC100PPSerialCommunication method</i> ), 142		set_init_shutter_status() ( <i>CryLasLaser method</i> ), 80
Sensor ( <i>class in hvl_ccb.utils.conversion.sensor</i> ), 176		set_jog_speed() ( <i>ILS2T method</i> ), 162
Sensor_error ( <i>PfeifferTPG.SensorStatus attribute</i> ), 144		set_lower_limits() ( <i>PSI9000 method</i> ), 87
Sensor_off ( <i>PfeifferTPG.SensorStatus attribute</i> ), 144		set_max_acceleration() ( <i>ILS2T method</i> ), 162
sensor_status ( <i>LuminoxMeasurementType attribute</i> ), 169		set_max_deceleration() ( <i>ILS2T method</i> ), 162
SERIAL ( <i>LaudaProRp245eConfig.ExtControlModeEnum attribute</i> ), 122		set_max_rpm() ( <i>ILS2T method</i> ), 162
serial_number ( <i>LuminoxMeasurementType attribute</i> ), 169		set_measuring_options() ( <i>MBW973 method</i> ), 126
SerialCommunication ( <i>class in hvl_ccb.comm.serial</i> ), 24		set_message_board() ( <i>BaseCube method</i> ), 38
SerialCommunicationBytesize ( <i>class in hvl_ccb.comm.serial</i> ), 25		set_motor_configuration() ( <i>NewportSMC100PP method</i> ), 136
SerialCommunicationConfig ( <i>class in hvl_ccb.comm.serial</i> ), 26		set_negative_software_limit() ( <i>NewportSMC100PP method</i> ), 136
SerialCommunicationIOError, 27		set_number_of_recordings() ( <i>HeinzingerDI method</i> ), 106
SerialCommunicationParity ( <i>class in hvl_ccb.comm.serial</i> ), 27		set_output() ( <i>PSI9000 method</i> ), 87
SerialCommunicationStopbits ( <i>class in hvl_ccb.comm.serial</i> ), 27		set_positive_software_limit() ( <i>NewportSMC100PP method</i> ), 136
Server ( <i>class in hvl_ccb.comm.opc</i> ), 24		set_pulse_energy() ( <i>CryLasLaser method</i> ), 80
		set_pump_level() ( <i>LaudaProRp245e method</i> ), 119
		set_ramp_iterations() ( <i>LaudaProRp245e method</i> ), 119
		set_ramp_program() ( <i>LaudaProRp245e method</i> ), 119
		set_ramp_segment() ( <i>LaudaProRp245e method</i> ), 119
		set_ramp_type() ( <i>ILS2T method</i> ), 162
		set_reference_point() ( <i>RTO1024 method</i> ), 156
		set_register() ( <i>FuGProbusV method</i> ), 97
		set_repetition_rate() ( <i>CryLasLaser method</i> ), 80

**set\_repetitions()** (*RTO1024 method*), 156  
**set\_status\_board()** (*BaseCube method*), 38  
**set\_system\_lock()** (*PSI9000 method*), 87  
**set\_temp\_set\_point()** (*LaudaProRp245e method*), 120  
**set\_transmission()** (*CryLasAttenuator method*), 76  
**set\_trigger\_level()** (*RTO1024 method*), 156  
**set\_trigger\_mode()** (*RTO1024 method*), 157  
**set\_trigger\_source()** (*RTO1024 method*), 157  
**set\_upper\_limits()** (*PSI9000 method*), 87  
**set\_voltage()** (*HeinzingerDI method*), 106  
**set\_voltage()** (*HeinzingerPNC method*), 108  
**set\_voltage\_current()** (*PSI9000 method*), 87  
**SetAttr** (*class in hvl\_ccb.utils.conversion.utils*), 177  
**SETCURRENT** (*FuGProbusVRegisterGroups attribute*), 100  
**setvalue** (*FuGProbusVSetRegisters property*), 100  
**SETVOLTAGE** (*FuGProbusVRegisterGroups attribute*), 100  
**SEVEN\_HUNDRED\_FIFTY\_VOLT** (*ACVoltageRange attribute*), 55  
**SEVENBITS** (*SerialCommunicationBytesize attribute*), 26  
**SHORT\_CIRCUIT** (*NewportSMC100PP.MotorErrors attribute*), 131  
**shunt** (*LEM4000S attribute*), 175  
**SHUTDOWN\_CURRENT\_LIMIT** (*PSI9000 attribute*), 86  
**SHUTDOWN\_VOLTAGE\_LIMIT** (*PSI9000 attribute*), 86  
**ShutterStatus** (*CryLasLaser attribute*), 79  
**ShutterStatus** (*CryLasLaserConfig attribute*), 82  
**SingleCommDevice** (*class in hvl\_ccb.dev.base*), 74  
**SIXBITS** (*SerialCommunicationBytesize attribute*), 26  
**SIXTEEN** (*HeinzingerConfig.RecordingsEnum attribute*), 104  
**SL** (*NewportConfigCommands attribute*), 130  
**SLOW\_FILTER** (*FilterRange attribute*), 56  
**SN** (*FuGReadbackChannels attribute*), 101  
**SOFTWARE\_INTERNAL\_SIXTY** (*CryLasLaser.RepetitionRates attribute*), 79  
**SOFTWARE\_INTERNAL\_TEN** (*CryLasLaser.RepetitionRates attribute*), 79  
**SOFTWARE\_INTERNAL\_TWENTY** (*CryLasLaser.RepetitionRates attribute*), 79  
**software\_revision** (*LuminoxMeasurementType attribute*), 169  
**SPACE** (*SerialCommunicationParity attribute*), 27  
**SPECIALRAMPUPWARDS** (*FuGRampModes attribute*), 101  
**spoll()** (*VisaCommunication method*), 33  
**spoll\_handler()** (*VisaDevice method*), 173  
**SR** (*NewportConfigCommands attribute*), 130  
**srq\_mask** (*FuGProbusVConfigRegisters property*), 97  
**srq\_status** (*FuGProbusVConfigRegisters property*), 97  
**stage\_configuration** (*NewportSMC100PPConfig attribute*), 139  
**START** (*LaudaProRp245eCommand attribute*), 121  
**start()** (*BaseCube method*), 38  
**start()** (*CryLasAttenuator method*), 76  
**start()** (*CryLasLaser method*), 81  
**start()** (*Device method*), 72  
**start()** (*DeviceSequenceMixin method*), 73  
**start()** (*ExperimentManager method*), 184  
**start()** (*Fluke8845a method*), 50  
**start()** (*FuG method*), 92  
**start()** (*FuGProbusIV method*), 96  
**start()** (*HeinzingerDI method*), 107  
**start()** (*HeinzingerPNC method*), 108  
**start()** (*ILS2T method*), 162  
**start()** (*LabJack method*), 117  
**start()** (*LaudaProRp245e method*), 120  
**start()** (*Luminox method*), 167  
**start()** (*MBW973 method*), 126  
**start()** (*NewportSMC100PP method*), 136  
**start()** (*PfeifferTPG method*), 147  
**start()** (*PSI9000 method*), 88  
**start()** (*RTO1024 method*), 157  
**start()** (*SingleCommDevice method*), 75  
**start()** (*Technix method*), 67  
**start()** (*VisaDevice method*), 173  
**start\_control()** (*MBW973 method*), 126  
**start\_polling()** (*Poller method*), 171  
**start\_ramp()** (*LaudaProRp245e method*), 120  
**STARTING** (*ExperimentStatus attribute*), 185  
**States** (*NewportSMC100PP attribute*), 131  
**status** (*BaseCube property*), 38  
**status** (*ExperimentManager property*), 184  
**status** (*FuGProbusVConfigRegisters property*), 98  
**status** (*FuGProbusVDORegisters property*), 99  
**status** (*Technix property*), 67  
**STATUSBYTE** (*FuGReadbackChannels attribute*), 101  
**STOP** (*LaudaProRp245eCommand attribute*), 121  
**stop()** (*BaseCube method*), 39  
**stop()** (*CryLasLaser method*), 81  
**stop()** (*Device method*), 72  
**stop()** (*DeviceSequenceMixin method*), 73  
**stop()** (*ExperimentManager method*), 184  
**stop()** (*Fluke8845a method*), 50  
**stop()** (*FuGProbusIV method*), 96  
**stop()** (*HeinzingerDI method*), 107  
**stop()** (*ILS2T method*), 162  
**stop()** (*LabJack method*), 117  
**stop()** (*LaudaProRp245e method*), 120  
**stop()** (*Luminox method*), 167  
**stop()** (*MBW973 method*), 126  
**stop()** (*NewportSMC100PP method*), 137  
**stop()** (*PfeifferTPG method*), 147  
**stop()** (*PSI9000 method*), 88  
**stop()** (*RTO1024 method*), 157  
**stop()** (*SingleCommDevice method*), 75  
**stop()** (*Technix method*), 67

**stop()** (*VisaDevice* method), 173  
**stop\_acquisition()** (*RTO1024* method), 157  
**stop\_motion()** (*NewportSMC100PP* method), 137  
**stop\_polling()** (*Poller* method), 171  
**stop\_ramp()** (*LaudaProRp245e* method), 120  
**STOP\_SAFETY\_STATUSES** (in module *hvl\_ccb.dev(cube.constants)*), 44  
**stopbits** (*CryLasAttenuatorSerialCommunicationConfig* attribute), 78  
**stopbits** (*CryLasLaserSerialCommunicationConfig* attribute), 85  
**stopbits** (*FuGSerialCommunicationConfig* attribute), 102  
**stopbits** (*HeinzingerSerialCommunicationConfig* attribute), 109  
**stopbits** (*LuminoxSerialCommunicationConfig* attribute), 171  
**stopbits** (*MBW973SerialCommunicationConfig* attribute), 129  
**stopbits** (*NewportSMC100PPSerialCommunicationConfig* attribute), 143  
**stopbits** (*PfeifferTPGSerialCommunicationConfig* attribute), 149  
**Stopbits** (*SerialCommunicationConfig* attribute), 26  
**stopbits** (*SerialCommunicationConfig* attribute), 27  
**streaming** (*LuminoxOutputMode* attribute), 169  
**StrEnumBase** (class in *hvl\_ccb.utils.enum*), 179  
**sub\_handler** (*BaseCubeOpcUaCommunicationConfig* attribute), 41  
**sub\_handler** (*OpcUaCommunicationConfig* attribute), 23  
**SyncCommunicationProtocol** (class in *hvl\_ccb.comm.base*), 15  
**SyncCommunicationProtocolConfig** (class in *hvl\_ccb.comm.base*), 15

**T**

**T** (*LabJack.ThermocoupleType* attribute), 113  
**t13\_socket\_1** (*BaseCube* attribute), 39  
**t13\_socket\_2** (*BaseCube* attribute), 39  
**t13\_socket\_3** (*BaseCube* attribute), 39  
**T1MS** (*FuGMonitorModes* attribute), 95  
**T200MS** (*FuGMonitorModes* attribute), 95  
**T20MS** (*FuGMonitorModes* attribute), 95  
**T256US** (*FuGMonitorModes* attribute), 95  
**T4** (*LabJack.DeviceType* attribute), 112  
**T4** (*LJMCommunicationConfig.DeviceType* attribute), 17  
**T40MS** (*FuGMonitorModes* attribute), 95  
**T4MS** (*FuGMonitorModes* attribute), 95  
**T560** (class in *hvl\_ccb.dev.highland\_t560.device*), 60  
**T560Communication** (class in *hvl\_ccb.dev.highland\_t560.base*), 58  
**T560CommunicationConfig** (class in *hvl\_ccb.dev.highland\_t560.base*), 58  
**T560Config** (class in *hvl\_ccb.dev.highland\_t560.device*), 61  
**T560Error**, 59  
**T7** (*LabJack.DeviceType* attribute), 112  
**T7** (*LJMCommunicationConfig.DeviceType* attribute), 17  
**T7\_PRO** (*LabJack.DeviceType* attribute), 112  
**T7\_PRO** (*LJMCommunicationConfig.DeviceType* attribute), 17  
**T800MS** (*FuGMonitorModes* attribute), 95  
**T80MS** (*FuGMonitorModes* attribute), 95  
**target\_pulse\_energy** (*CryLasLaser* property), 81  
**Tcp** (class in *hvl\_ccb.comm.tcp*), 28  
**TCP** (*LJMCommunicationConfig.ConnectionType* attribute), 17  
**TcpCommunicationConfig** (class in *hvl\_ccb.comm.tcp*), 28  
**TCPIP\_INSTR** (*VisaCommunicationConfig.InterfaceType* attribute), 33  
**TCPIP\_SOCKET** (*VisaCommunicationConfig.InterfaceType* attribute), 33  
**TEC1** (*CryLasLaser.AnswersStatus* attribute), 78  
**TEC2** (*CryLasLaser.AnswersStatus* attribute), 78  
**Technix** (class in *hvl\_ccb.dev.technix.device*), 66  
**TechnixConfig** (class in *hvl\_ccb.dev.technix.device*), 67  
**TechnixError**, 63  
**TechnixFaultError**, 63  
**TechnixSerialCommunication** (class in *hvl\_ccb.dev.technix.base*), 63  
**TechnixSerialCommunicationConfig** (class in *hvl\_ccb.dev.technix.base*), 63  
**TechnixTelnetCommunication** (class in *hvl\_ccb.dev.technix.base*), 64  
**TechnixTelnetCommunicationConfig** (class in *hvl\_ccb.dev.technix.base*), 64  
**TelnetCommunication** (class in *hvl\_ccb.comm.telnet*), 30  
**TelnetCommunicationConfig** (class in *hvl\_ccb.comm.telnet*), 31  
**TelnetError**, 32  
**TEMP** (*ILS2TRegAddr* attribute), 165  
**TEMP\_SET\_POINT** (*LaudaProRp245eCommand* attribute), 121  
**temp\_set\_point\_init** (*LaudaProRp245eConfig* attribute), 123  
**Temperature** (class in *hvl\_ccb.utils.conversion.unit*), 177  
**temperature\_sensor** (*LuminoxMeasurementType* attribute), 169  
**temperature\_unit** (*LMT70A* attribute), 176  
**TEN** (*LabJack.AInRange* attribute), 111  
**TEN** (*LabJack.CalMicroAmpere* attribute), 111  
**TEN\_AMPERE** (*ACCurrrentRange* attribute), 54  
**TEN\_AMPERE** (*DCCurrentRange* attribute), 55  
**TEN\_MHZ** (*LabJack.ClockFrequency* attribute), 112

TEN\_MILLI\_AMPERE (*ACCurrerntRange attribute*), 54  
 TEN\_MILLI\_AMPERE (*DCCurrentRange attribute*), 55  
 TEN\_MILLI\_SECOND (*ApertureRange attribute*), 55  
 TEN\_MILLION\_OHM (*ResistanceRange attribute*), 56  
 TEN\_THOUSAND\_OHM (*ResistanceRange attribute*), 56  
 TEN\_VOLT (*ACVoltageRange attribute*), 55  
 TEN\_VOLT (*DCVoltageRange attribute*), 56  
**terminator** (*AsyncCommunicationProtocolConfig attribute*), 14  
**terminator** (*CryLasAttenuatorSerialCommunicationConfig attribute*), 78  
**terminator** (*CryLasLaserSerialCommunicationConfig attribute*), 85  
**terminator** (*Fluke8845aTelnetCommunicationConfig attribute*), 52  
**TERMINATOR** (*FuGProbusIVCommands attribute*), 96  
**terminator** (*FuGProbusVConfigRegisters property*), 98  
**terminator** (*FuGSerialCommunicationConfig attribute*), 103  
**terminator** (*HeinzingerSerialCommunicationConfig attribute*), 110  
**terminator** (*LaudaProRp245eTcpCommunicationConfig attribute*), 125  
**terminator** (*LuminoxSerialCommunicationConfig attribute*), 171  
**terminator** (*MBW973SerialCommunicationConfig attribute*), 129  
**terminator** (*NewportSMC100PPSerialCommunicationConfig attribute*), 143  
**terminator** (*PfeifferTPGSerialCommunicationConfig attribute*), 150  
**terminator** (*T560CommunicationConfig attribute*), 59  
**terminator\_str()** (*SerialCommunicationConfig method*), 27  
 THIRTY\_TWO\_BIT (*LabJack.BitLimit attribute*), 111  
 THREE\_AMPERE (*ACCurrerntRange attribute*), 54  
 THREE\_AMPERE (*DCCurrentRange attribute*), 55  
**timeout** (*CryLasAttenuatorSerialCommunicationConfig attribute*), 78  
**timeout** (*CryLasLaserSerialCommunicationConfig attribute*), 85  
**timeout** (*FuGSerialCommunicationConfig attribute*), 103  
**timeout** (*HeinzingerSerialCommunicationConfig attribute*), 110  
**timeout** (*LuminoxSerialCommunicationConfig attribute*), 171  
**timeout** (*MBW973SerialCommunicationConfig attribute*), 129  
**timeout** (*NewportSMC100PPSerialCommunicationConfig attribute*), 143  
**timeout** (*PfeifferTPGSerialCommunicationConfig attribute*), 150  
**timeout** (*SerialCommunicationConfig attribute*), 27  
**timeout** (*TelnetCommunicationConfig attribute*), 31  
**timeout** (*VisaCommunicationConfig attribute*), 35  
**timeout\_interval** (*BaseCubeConfiguration attribute*), 40  
**timeout\_status\_change** (*BaseCubeConfiguration attribute*), 40  
**timeout\_test\_parameters** (*PICubeConfiguration attribute*), 47  
**TORR** (*Pressure attribute*), 177  
 TPG25xA (*PfeifferTPGConfig.Model attribute*), 147  
 TPGx6x (*PfeifferTPGConfig.Model attribute*), 147  
**TPR** (*PfeifferTPG.SensorTypes attribute*), 145  
**transmission** (*CryLasAttenuator property*), 76  
**trigger()** (*Fluke8845a method*), 50  
**trigger\_delay** (*Fluke8845a property*), 50  
**trigger\_level** (*T560 property*), 61  
**trigger\_mode** (*T560 property*), 61  
**trigger\_source** (*Fluke8845a property*), 50  
**TriggerMode** (*class in hvl\_ccb.dev.hIGHLAND\_t560.base*), 59  
**TriggerSource** (*class in hvl\_ccb.dev.fluke884x.constants*), 54  
**TWELVE\_HUNDRED\_FIFTY\_KHZ** (*LabJack.ClockFrequency attribute*), 112  
**TWENTY\_FIVE\_HUNDRED\_KHZ** (*LabJack.ClockFrequency attribute*), 112  
**TWENTY\_MHZ** (*LabJack.ClockFrequency attribute*), 112  
**TWO** (*HeinzingerConfig.RecordingsEnum attribute*), 104  
**TWO** (*SerialCommunicationStopbits attribute*), 28  
**TWO\_HUNDRED** (*LabJack.CalMicroAmpere attribute*), 111  
**TWO\_WIRE\_RESISTANCE** (*MeasurementFunction attribute*), 53  
**two\_wire\_resistance\_range** (*Fluke8845a attribute*), 51

## U

**Underrange** (*PfeifferTPG.SensorStatus attribute*), 145  
**Unit** (*class in hvl\_ccb.utils.conversion.unit*), 177  
**unit** (*ILS2TModbusTcpCommunicationConfig attribute*), 165  
**unit** (*ModbusTcpCommunicationConfig attribute*), 20  
**unit** (*PfeifferTPG property*), 147  
**unit()** (*RangeEnum class method*), 179  
**unit\_current** (*HeinzingerPNC property*), 108  
**unit\_voltage** (*HeinzingerPNC property*), 108  
**UNKNOWN** (*HeinzingerDI.OutputStatus attribute*), 105  
**UNKNOWN** (*HeinzingerPNC.UnitCurrent attribute*), 107  
**UNKNOWN** (*HeinzingerPNC.UnitVoltage attribute*), 107  
**UNREADY\_INACTIVE** (*CryLasLaser.LaserStatus attribute*), 79  
**update\_laser\_status()** (*CryLasLaser method*), 81  
**update\_parameter** (*OpcUaCommunicationConfig attribute*), 23

**update\_repetition\_rate()** (*CryLasLaser method*), 81  
**update\_shutter\_status()** (*CryLasLaser method*), 81  
**update\_target\_pulse\_energy()** (*CryLasLaser method*), 81  
**UpdateEspStageInfo** (*NewportSMC100PPConfig.EspStageConfig attribute*), 138  
**UPPER\_TEMP** (*LaudaProRp245eCommand attribute*), 121  
**upper\_temp** (*LaudaProRp245eConfig attribute*), 123  
**USB** (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 122  
**USB** (*LJMCommunicationConfig.ConnectionType attribute*), 17  
**use\_external\_clock()** (*T560 method*), 61  
**user\_position\_offset** (*NewportSMC100PPConfig attribute*), 139  
**user\_steps()** (*ILS2T method*), 163

**V**

**V** (*HeinzingerPNC.UnitVoltage attribute*), 107  
**VA** (*NewportConfigCommands attribute*), 130  
**validate\_and\_resolve\_host()** (*in module hvl\_ccb.utils.validation*), 180  
**validate\_bool()** (*in module hvl\_ccb.utils.validation*), 180  
**validate\_number()** (*in module hvl\_ccb.utils.validation*), 180  
**validate\_pump\_level()** (*LaudaProRp245e method*), 120  
**validate\_tcp\_port()** (*in module hvl\_ccb.utils.validation*), 181  
**value** (*FuGProbusVMonitorRegisters property*), 99  
**value** (*LabJack.AInRange property*), 111  
**value\_raw** (*FuGProbusVMonitorRegisters property*), 99  
**ValueEnum** (*class in hvl\_ccb.utils.enum*), 179  
**VB** (*NewportConfigCommands attribute*), 130  
**velocity** (*NewportSMC100PPConfig attribute*), 139  
**visa\_backend** (*VisaCommunicationConfig attribute*), 35  
**VisaCommunication** (*class in hvl\_ccb.comm.visa*), 32  
**VisaCommunicationConfig** (*class in hvl\_ccb.comm.visa*), 33  
**VisaCommunicationConfig.InterfaceType** (*class in hvl\_ccb.comm.visa*), 33  
**VisaCommunicationError**, 35  
**VisaDevice** (*class in hvl\_ccb.dev.visa*), 172  
**VisaDeviceConfig** (*class in hvl\_ccb.dev.visa*), 173  
**VOLT** (*ILS2TRegAddr attribute*), 165  
**voltage** (*FuG property*), 92  
**VOLTAGE** (*FuGProbusIVCommands attribute*), 97  
**VOLTAGE** (*FuGReadbackChannels attribute*), 101  
**voltage** (*Technix property*), 67  
**VOLTAGE\_AC** (*MeasurementFunction attribute*), 53  
**voltage\_actual** (*PICube property*), 46  
**VOLTAGE\_DC** (*MeasurementFunction attribute*), 53  
**voltage\_filter** (*Fluke8845a attribute*), 51  
**voltage\_lower\_limit** (*PSI9000Config attribute*), 89  
**voltage\_max** (*PICube property*), 46  
**voltage\_monitor** (*FuG property*), 92  
**voltage\_primary** (*PICube property*), 46  
**voltage\_regulation** (*Technix property*), 67  
**voltage\_upper\_limit** (*PSI9000Config attribute*), 89

**W**

**WAIT\_AFTER\_WRITE** (*VisaCommunication attribute*), 32  
**wait\_for\_polling\_result()** (*Poller method*), 171  
**wait\_operation\_complete()** (*VisaDevice method*), 173  
**wait\_sec\_initialisation** (*PSI9000Config attribute*), 89  
**wait\_sec\_max\_disable** (*ILS2TConfig attribute*), 164  
**wait\_sec\_post\_absolute\_position** (*ILS2TConfig attribute*), 164  
**wait\_sec\_post\_activate** (*LuminoxConfig attribute*), 168  
**wait\_sec\_post\_cannot\_disable** (*ILS2TConfig attribute*), 164  
**wait\_sec\_post\_enable** (*ILS2TConfig attribute*), 164  
**wait\_sec\_post\_relative\_step** (*ILS2TConfig attribute*), 164  
**wait\_sec\_pre\_read\_or\_write** (*LaudaProRp245eTcpCommunicationConfig attribute*), 125  
**wait\_sec\_read\_text\_nonempty** (*AsyncCommunicationProtocolConfig attribute*), 14  
**wait\_sec\_read\_text\_nonempty** (*FuGSerialCommunicationConfig attribute*), 103  
**wait\_sec\_read\_text\_nonempty** (*HeinzingerSerialCommunicationConfig attribute*), 110  
**wait\_sec\_settings\_effect** (*PSI9000Config attribute*), 89  
**wait\_sec\_stop\_commands** (*FuGConfig attribute*), 93  
**wait\_sec\_stop\_commands** (*HeinzingerConfig attribute*), 104  
**wait\_sec\_system\_lock** (*PSI9000Config attribute*), 89  
**wait\_sec\_trials\_activate** (*LuminoxConfig attribute*), 168  
**wait\_timeout\_retry\_sec** (*OpcUaCommunicationConfig attribute*), 23  
**wait\_until\_motor\_initialized()** (*NewportSMC100PP method*), 137  
**wait\_until\_ready()** (*CryLasLaser method*), 81  
**WIFI** (*LJMCommunicationConfig.ConnectionType attribute*), 17  
**write()** (*AsyncCommunicationProtocol method*), 13  
**write()** (*BaseCube method*), 39  
**write()** (*MBW973 method*), 127  
**write()** (*OpcUaCommunication method*), 22

`write()` (*Tcp method*), 28  
`write()` (*VisaCommunication method*), 33  
`write_absolute_position()` (*ILS2T method*), 163  
`write_bytes()` (*AsyncCommunicationProtocol method*), 13  
`write_bytes()` (*SerialCommunication method*), 25  
`write_bytes()` (*TelnetCommunication method*), 30  
`write_command()` (*LaudaProRp245eTcpCommunication method*), 124  
`write_name()` (*LJMCommunication method*), 17  
`write_names()` (*LJMCommunication method*), 17  
`write_registers()` (*ModbusTcpCommunication method*), 20  
`write_relative_step()` (*ILS2T method*), 163  
`write_termination` (*VisaCommunicationConfig attribute*), 35  
`write_text()` (*AsyncCommunicationProtocol method*), 13  
`WRONG_ESP_STAGE` (*NewportSMC100PP.MotorErrors attribute*), 131

## X

`x_stat` (*FuGProbusVDIRegisters property*), 98  
`XOUTPUTS` (*FuGProbusIVCommands attribute*), 97

## Y

`YES` (*FuGDigitalVal attribute*), 93

## Z

`ZX` (*NewportConfigCommands attribute*), 130