

---

# HVL Common Code Base Documentation

*Release 0.6.1*

**Mikolaj Rybiński, David Graber, Henrik Menne, Alise Chachereau,**

**May 08, 2021**



# CONTENTS:

<b>1</b>	<b>HVL Common Code Base</b>	<b>1</b>
1.1	Features . . . . .	1
1.2	Documentation . . . . .	4
1.3	Credits . . . . .	4
<b>2</b>	<b>Installation</b>	<b>5</b>
2.1	Stable release . . . . .	5
2.2	From sources . . . . .	5
2.3	Additional system libraries . . . . .	6
<b>3</b>	<b>Usage</b>	<b>7</b>
<b>4</b>	<b>API Documentation</b>	<b>9</b>
4.1	hvl_ccb . . . . .	9
<b>5</b>	<b>Contributing</b>	<b>169</b>
5.1	Types of Contributions . . . . .	169
5.2	Get Started! . . . . .	170
5.3	Merge Request Guidelines . . . . .	171
5.4	Tips . . . . .	171
5.5	Deploying . . . . .	172
<b>6</b>	<b>Credits</b>	<b>173</b>
6.1	Maintainers . . . . .	173
6.2	Authors . . . . .	173
6.3	Contributors . . . . .	173
<b>7</b>	<b>History</b>	<b>175</b>
7.1	0.6.1 (2021-05-08) . . . . .	175
7.2	0.6.0 (2021-04-23) . . . . .	175
7.3	0.5.0 (2020-11-11) . . . . .	176
7.4	0.4.0 (2020-07-16) . . . . .	176
7.5	0.3.5 (2020-02-18) . . . . .	177
7.6	0.3.4 (2019-12-20) . . . . .	177
7.7	0.3.3 (2019-05-08) . . . . .	178
7.8	0.3.2 (2019-05-08) . . . . .	178
7.9	0.3.1 (2019-05-02) . . . . .	178
7.10	0.3 (2019-05-02) . . . . .	178
7.11	0.2.1 (2019-04-01) . . . . .	178
7.12	0.2.0 (2019-03-31) . . . . .	178
7.13	0.1.0 (2019-02-06) . . . . .	179

<b>8 Indices and tables</b>	<b>181</b>
<b>Python Module Index</b>	<b>183</b>
<b>Index</b>	<b>185</b>

---

**CHAPTER  
ONE**

---

## **HVL COMMON CODE BASE**

Python common code base to control devices high voltage research devices, in particular, as used in Christian Franck's High Voltage Lab (HVL), D-ITET, ETH.

- Free software: GNU General Public License v3
- Copyright (c) 2019-2021 ETH Zurich, SIS ID and HVL D-ITET

### **1.1 Features**

For managing multi-device experiments instantiate the `ExperimentManager` utility class.

#### **1.1.1 Devices**

The devices wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, various settings and options enumerations, as well as start/stop methods. Currently, wrappers to control the following devices are available:

Function/Type	Devices
Data acquisition and Digital IO	LabJack (T4, T7, T7-PRO)
Experiment control	HVL Supercube with and without Frequency Converter
Gas Analyser	MBW 973-SF6 gas dew point mirror analyzer Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges SST Luminox oxygen sensor
I2C host	TiePie (HS5, WS5)
Laser	CryLaS pulsed laser CryLaS laser attenuator
Oscilloscope	Rhode & Schwarz RTO 1024 TiePie (HS5, HS6, WS5)
Power supply	Elektro-Automatik PSI9000 FuG Elektronik Heinzinger PNC Technix capacitor charger
Stepper motor drive	Newport SMC100PP Schneider Electric ILS2T
Waveform generator	TiePie (HS5, WS5)

Each device uses at least one standardised communication protocol wrapper.

### 1.1.2 Communication protocols

In hvl\_ccb by “communication protocol” we mean different levels of communication standards, from the low level actual communication protocols like serial communication to application level interfaces like VISA TCP standard. There are also devices in hvl\_ccb that use dummy communication protocol concept; this is because these devices build on propriety vendor libraries that communicate with vendor devices, like in case of the TiePie devices.

The communication protocol wrappers in hvl\_ccb provide a standardised API with configuration dataclasses, as well as open/close, and read/write/query methods. Currently, wrappers to use the following communication protocols are available:

Communication protocol	Devices using
Modbus TCP	Schneider Electric ILS2T stepper motor drive
OPC UA	HVL Supercube with and without Frequency Converter
Serial	CryLaS pulsed laser and laser attenuator FuG Elektronik power supply (e.g. capacitor charger HCK) using the Probus V protocol Heinzinger PNC power supply using Heinzinger Digital Interface I/II SST Luminox oxygen sensor MBW 973-SF6 gas dew point mirror analyzer Newport SMC100PP single axis driver for 2-phase stepper motors Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges Technix capacitor charger
Telnet	Technix capacitor charger
VISA TCP	Elektro-Automatik PSI9000 DC power supply Rhode & Schwarz RTO 1024 oscilloscope
<i>proprietary</i>	LabJack (T4, T7, T7-PRO) devices, which communicate via LJM Library TiePie (HS5, HS6, WS5) oscilloscopes, generators and I2C hosts, which communicate via LibTiePie SDK

## 1.2 Documentation

Note: if you're planning to contribute to the hvl\_ccb project do read beforehand the **Contributing** section in the HVL CCB documentation.

Do either:

- read [HVL CCB documentation at RTD](#),

or

- build and read HVL CCB documentation locally; install first documentation build requirements:

```
$ pip install docs/requirements.txt
```

and then either on Windows in Git BASH run:

```
$ ./make.sh docs
```

or from any other shell with GNU Make installed run:

```
$ make docs
```

The target index HTML ("docs/\_build/html/index.html") should open automatically in your Web browser.

## 1.3 Credits

This package was created with [Cookiecutter](#) and the [audreyr/cookiecutter-pypackage](#) project template.

---

CHAPTER  
TWO

---

## INSTALLATION

### 2.1 Stable release

To install HVL Common Code Base, run this command in your terminal:

```
$ pip install hvl_ccb
```

To install HVL Common Code Base with optional Python libraries that require additional external libraries installations like tiepie or labjack, specify on installation extra features you want by running e.g. this command in your terminal:

```
$ pip install hvl_ccb[tiepie,labjack]
```

This is the preferred method to install HVL Common Code Base, as it will always install the most recent stable release.

If you don't have `pip` installed, this [Python installation guide](#) can guide you through the process.

### 2.2 From sources

The sources for HVL Common Code Base can be downloaded from the [GitLab](#) repo.

You can either clone the repository:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

Or download the [tarball](#):

```
$ curl -OL https://gitlab.com/ethz_hvl/hvl_ccb/-/archive/master/hvl_ccb.tar.gz
```

Once you have a copy of the source, you can install it with:

```
$ pip install .
```

## 2.3 Additional system libraries

Please note that for some of the dependencies, like e.g. for `labjack-ljm` or `python-libtiepie`, you may need to separately install additional system libraries.

For [LJM Library](#) please make sure that you have installed version 2019\_02\_14 or higher.

---

**CHAPTER  
THREE**

---

**USAGE**

To use HVL Common Code Base in a project:

```
import hvl_ccb
```



## API DOCUMENTATION


### 4.1 hvl\_ccb

#### 4.1.1 Subpackages

[hvl\\_ccb.comm](#)

**Submodules**

[hvl\\_ccb.comm.base](#)

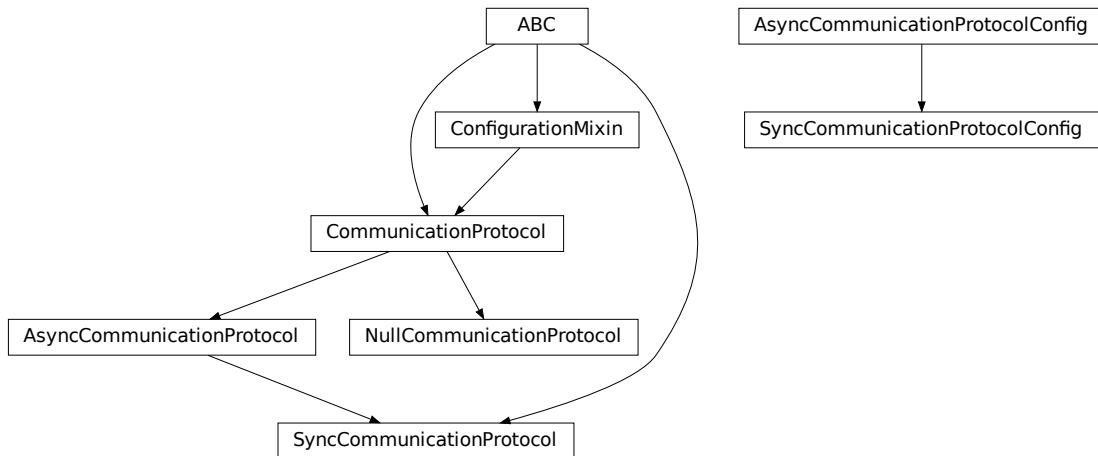

Module with base classes for communication protocols.

**class** `AsyncCommunicationProtocol` (`config`)  
Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Abstract base class for asynchronous communication protocols

**static config\_cls()** → Type[`hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`]  
Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

### `read()` → str

Read a single line of text as *str* from the communication.

**Returns** text as *str* including the terminator, which can also be empty “”

### `read_all(n_attempts_max: Optional[int] = None, attempt_interval_sec: Optional[Union[int, float]] = None)` → Optional[str]

Read all lines of text from the connection till nothing is left to read.

#### Parameters

- **n\_attempts\_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt\_interval\_sec** – time between the reading attempts

**Returns** A multi-line *str* including the terminator internally

### `abstract read_bytes()` → bytes

Read a single line as *bytes* from the communication.

This method uses *self.access\_lock* to ensure thread-safety.

**Returns** a single line as *bytes* containing the terminator, which can also be empty b””

### `read_nonempty(n_attempts_max: Optional[int] = None, attempt_interval_sec: Optional[Union[int, float]] = None)` → Optional[str]

Try to read a non-empty single line of text as *str* from the communication. If the host does not reply or reply with white space only, it will return None.

**Returns** a non-empty text as a *str* or *None* in case of an empty string

#### Parameters

- **n\_attempts\_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt\_interval\_sec** – time between the reading attempts

### `read_text()` → str

Read one line of text from the serial port. The input buffer may hold additional data afterwards, since only one line is read.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

**Returns** String read from the serial port; “” if there was nothing to read.

**Raises** `SerialCommunicationIOError` – when communication port is not opened

### `read_text_nonempty(n_attempts_max: Optional[int] = None, attempt_interval_sec: Optional[Union[int, float]] = None)` → Optional[str]

Reads from the serial port, until a non-empty line is found, or the number of attempts is exceeded.

NOTE: backward-compatibility proxy for *read\_text* method; to be removed in v1.0

Attention: in contrast to *read\_text*, the returned answer will be stripped of a whitespace newline terminator at the end, if such terminator is set in the initial configuration (default).

#### Parameters

- **n\_attempts\_max** – maximum number of read attempts
- **attempt\_interval\_sec** – time between the reading attempts

**Returns** String read from the serial port; “” if number of attempts is exceeded or serial port is not opened.

### `write(text: str)`

Write text as *str* to the communication.

**Parameters** `text` – test as a `str` to be written

**abstract** `write_bytes` (`data: bytes`) → `int`  
Write data as `bytes` to the communication.  
This method uses `self.access_lock` to ensure thread-safety.

**Parameters** `data` – data as `bytes`-string to be written

**Returns** number of bytes written

`write_text` (`text: str`)  
Write text to the serial port. The text is encoded and terminated by the configured terminator.  
NOTE: backward-compatibility proxy for `read` method; to be removed in v1.0

**Parameters** `text` – Text to send to the port.

**Raises** `SerialCommunicationIOError` – when communication port is not opened

```
class AsyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding_error_handling: str = 'replace', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10)
```

Bases: `object`

Base configuration data class for asynchronous communication protocols

`clean_values()`

`default_n_attempts_read_text_nonempty: int = 10`  
default number of attempts to read a non-empty text

`encoding: str = 'utf-8'`  
Standard encoding of the connection. Typically this is `utf-8`, but can also be `latin-1` or something from here: <https://docs.python.org/3/library/codecs.html#standard-encodings>

`encoding_error_handling: str = 'replace'`  
Encoding error handling scheme as defined here: <https://docs.python.org/3/library/codecs.html#error-handlers> By default replacing invalid characters with “`UFFFD`” REPLACEMENT CHARACTER on decoding and with “`?`” on decoding.

`force_value` (`fieldname, value`)  
Forces a value to a dataclass field despite the class being frozen.  
NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- `fieldname` – name of the field
- `value` – value to assign

`is_configdataclass = True`

`classmethod keys()` → `Sequence[str]`  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

`classmethod optional_defaults()` → `Dict[str, object]`  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**terminator: bytes = b'\r\n'**

The terminator character. Typically this is `b'\r\n'` or `b'\n'`, but can also be `b'\r'` or other combinations. This defines the end of a single line.

**wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5**

time to wait between attempts of reading a non-empty text

**class CommunicationProtocol(config)**

Bases: `hvl_ccb.configuration.ConfigurationMixin`, `abc.ABC`

Communication protocol abstract base class.

Specifies the methods to implement for communication protocol, as well as implements some default settings and checks.

**access\_lock**

Access lock to use with context manager when accessing the communication protocol (thread safety)

**abstract close()**

Close the communication protocol

**abstract open()**

Open communication protocol

**class NullCommunicationProtocol(config)**

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Communication protocol that does nothing.

**close()** → None

Void close function.

**static config\_cls()** → Type[`hvl_ccb.configuration.EmptyConfig`]

Empty configuration

**Returns** EmptyConfig

**open()** → None

Void open function.

**class SyncCommunicationProtocol(config)**

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocol`, `abc.ABC`

Abstract base class for synchronous communication protocols with `query()`

**static config\_cls()** → Type[`hvl_ccb.comm.base.SyncCommunicationProtocolConfig`]

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query(command: str)** → Optional[str]


Send a command to the interface and handle the status message. Eventually raises an exception.

**Parameters** `command` – Command to send

**Returns** Answer from the interface, which can be None instead of an empty reply

```
class SyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str =
    'utf-8', encoding_error_handling: str = 'replace',
    wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int =
    10)
Bases: hvl_ccb.comm.base.AsyncCommunicationProtocolConfig
```

## hvl\_ccb.comm.labjack\_ljm

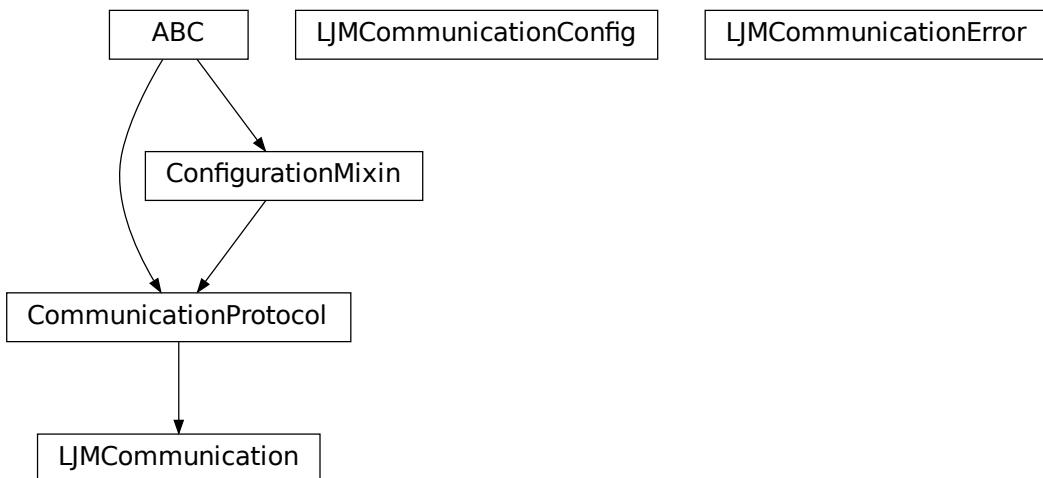

Communication protocol for LabJack using the LJM Library. Originally developed and tested for LabJack T7-PRO.

Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

```
class LJMCommunication(configuration)
Bases: hvl_ccb.comm.base.CommunicationProtocol
```

Communication protocol implementing the LabJack LJM Library Python wrapper.

**close()** → None  
Close the communication port.

**static config\_cls()**  
Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property is\_open**  
Flag indicating if the communication port is open.

**Returns** `True` if the port is open, otherwise `False`

**open()** → None  
Open the communication port.

```
read_name(*names: str, return_num_type: Type[numbers.Real] = <class 'float'>) → Union[numbers.Real, Sequence[numbers.Real]]
```

Read one or more input numeric values by name.

#### Parameters

- **names** – one or more names to read out from the LabJack
- **return\_num\_type** – optional numeric type specification for return values; by default *float*.

**Returns** answer of the LabJack, either single number or multiple numbers in a sequence, respectively, when one or multiple names to read were given

**Raises** **TypeError** – if read value of type not compatible with *return\_num\_type*

```
write_name(name: str, value: numbers.Real) → None
```

Write one value to a named output.

#### Parameters

- **name** – String or with name of LabJack IO
- **value** – is the value to write to the named IO port

```
write_names(name_value_dict: Dict[str, numbers.Real]) → None
```

Write more than one value at once to named outputs.

**Parameters** **name\_value\_dict** – is a dictionary with string names of LabJack IO as keys and corresponding numeric values

```
class LJMCommunicationConfig(device_type: Union[str, hvl_ccb._dev.labjack.DeviceType]
                                = 'ANY', connection_type: Union[str, hvl_ccb.comm.labjack_ljm.LJMCCommunicationConfig.ConnectionType]
                                = 'ANY', identifier: str = 'ANY')
```

Bases: *object*

Configuration dataclass for *LJMCommunication*.

```
class ConnectionType(value=<object object>, names=None, module=None, type=None,
                      start=1, boundary=None)
```

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

LabJack connection type.

**ANY** = 1

**ETHERNET** = 4

**TCP** = 3

**USB** = 2

**WIFI** = 5

```
class DeviceType(value=<object object>, names=None, module=None, type=None, start=1,
                    boundary=None)
```

Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

LabJack device types.

Can be also looked up by ambiguous Product ID (*p\_id*) or by instance name: `python LabJackDeviceType(4)` is *LabJackDeviceType('T4')* `

**ANY** = 1

**T4** = 2

```

T7 = 3
T7_PRO = 4

classmethod get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
List[hvl_ccb._dev.labjack.DeviceType]]
Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

Parameters p_id – Product ID of a LabJack device
Returns Instance or list of instances of LabJackDeviceType
Raises ValueError – when Product ID is unknown

clean_values() → None
Performs value checks on device_type and connection_type.

connection_type: Union[str, hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType]
Can be either string or of enum ConnectionType.

device_type: Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY'
Can be either string 'ANY', 'T7_PRO', 'T7', 'T4', or of enum DeviceType.

force_value(fieldname, value)
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define post_force_value method with same signature as this method to do extra processing
after value has been forced on fieldname.

Parameters

- fieldname – name of the field
- value – value to assign

identifier: str = 'ANY'
The identifier specifies information for the connection to be used. This can be an IP address, serial number,
or device name. See the LabJack docs ( https://labjack.com/support/software/api/ljm/function-reference/ljmopens/identifier-parameter) for more information.

is_configdataclass = True

classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
instantiation.


Returns a list of strings containing all required keys.

exception LJMCommunicationError
Bases: Exception

Errors coming from LJMCommunication.

```

## `hvl_ccb.comm.modbus_tcp`

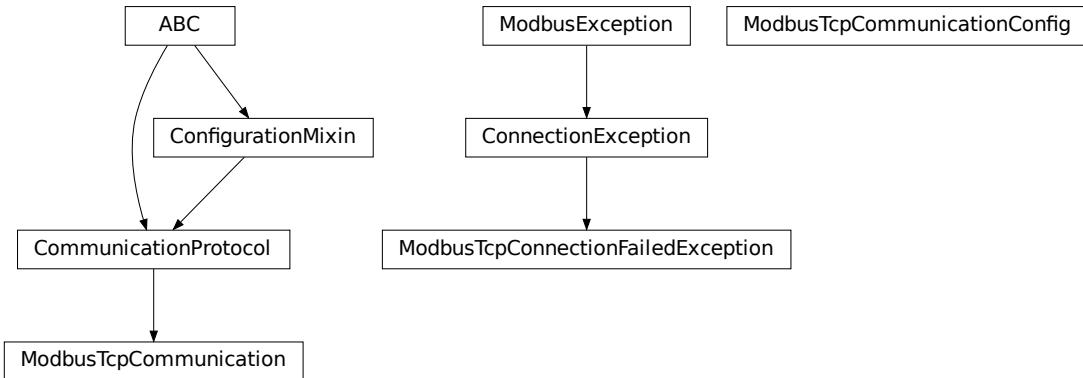

Communication protocol for modbus TCP ports. Makes use of the `pymodbus` library.

**class** `ModbusTcpCommunication`(*configuration*)  
*Bases:* `hvl_ccb.comm.base.CommunicationProtocol`

Implements the Communication Protocol for modbus TCP.

**close()**

Close the Modbus TCP connection.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**open()** → None

Open the Modbus TCP connection.

**Raises** `ModbusTcpConnectionFailedException` – if the connection fails.

**read\_holding\_registers**(*address: int, count: int*) → List[int]

Read specified number of register starting with given address and return the values from each register.

**Parameters**

- **address** – address of the first register
- **count** – count of registers to read

**Returns** list of *int* values

**read\_input\_registers**(*address: int, count: int*) → List[int]

Read specified number of register starting with given address and return the values from each register in a list.

**Parameters**

- **address** – address of the first register
- **count** – count of registers to read

**Returns** list of *int* values

---

**write\_registers** (*address*: int, *values*: Union[List[int], int])

Write values from the specified address forward.

#### Parameters

- **address** – address of the first register
- **values** – list with all values

**class ModbusTcpCommunicationConfig** (*host*: str, *unit*: int, *port*: int = 502)

Bases: object

Configuration dataclass for *ModbusTcpCommunication*.

**clean\_values** ()

**force\_value** (*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**host**: str

Host is the IP address of the connected device.

**is\_configdataclass** = True

**classmethod keys** () → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults** () → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port**: int = 502

TCP port

**classmethod required\_keys** () → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**unit**: int


Unit number to be used when connecting with Modbus/TCP. Typically this is used when connecting to a relay having Modbus/RTU-connected devices.

**exception ModbusTcpConnectionFailedException** (*string*=")

Bases: pymodbus.exceptions.ConnectionException

Exception raised when the connection failed.

## hvl\_ccb.comm.opc

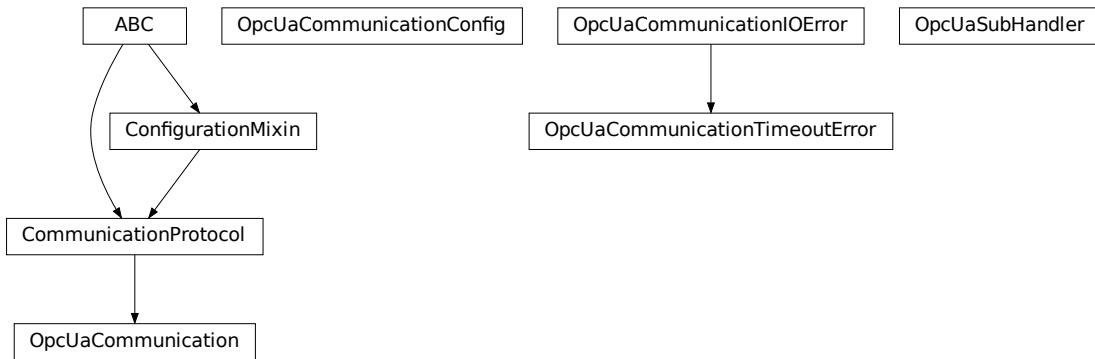

Communication protocol implementing an OPC UA connection. This protocol is used to interface with the “Super-cube” PLC from Siemens.

**class OpcUaCommunication(config)**  
Bases: [hvl\\_ccb.comm.base.CommunicationProtocol](#)

Communication protocol implementing an OPC UA connection. Makes use of the package python-opcua.

**close() → None**

Close the connection to the OPC UA server.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**init\_monitored\_nodes(node\_id: Union[object, Iterable], ns\_index: int) → None**

Initialize monitored nodes.

### Parameters

- **node\_id** – one or more strings of node IDs; node IDs are always casted via `str()` method here, hence do not have to be strictly string objects.
- **ns\_index** – the namespace index the nodes belong to.

**Raises** `OpcUaCommunicationIOError` – when protocol was not opened or can't communicate with a OPC UA server

**property is\_open**

Flag indicating if the communication port is open.

**Returns** `True` if the port is open, otherwise `False`

**open() → None**

Open the communication to the OPC UA server.

**Raises** `OpcUaCommunicationIOError` – when communication port cannot be opened.

**read(node\_id, ns\_index)**

Read a value from a node with id and namespace index.

### Parameters

- **node\_id** – the ID of the node to read the value from
- **ns\_index** – the namespace index of the node

**Returns** the value of the node object.

**Raises** *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

**write**(*node\_id*, *ns\_index*, *value*) → None

Write a value to a node with name *name*.

#### Parameters

- **node\_id** – the id of the node to write the value to.
- **ns\_index** – the namespace index of the node.
- **value** – the value to write.

**Raises** *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

```
class OpcUaCommunicationConfig(host: str, endpoint_name: str, port: int = 4840,
                                sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
                                <hvl_ccb.comm.opc.OpcUaSubHandler object>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)
```

Bases: *object*

Configuration dataclass for OPC UA Communciation.

**clean\_values**()

**endpoint\_name**: str

Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**host**: str

Hostname or IP-Address of the OPC UA server.

**is\_configdataclass** = True

**classmethod keys**() → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**max\_timeout\_retry\_nr**: int = 5

Maximal number of call re-tries on underlying OPC UA client timeout error

**classmethod optional\_defaults**() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```
port: int = 4840
Port of the OPC UA server to connect to.

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.comm.opc.OpcUaSubHandler object
object to use for handling subscriptions.

update_period: int = 500
Update period for generating datachange events in OPC UA [milli seconds]

wait_timeout_retry_sec: Union[int, float] = 1
Wait time between re-trying calls on underlying OPC UA client timeout error

exception OpcUaCommunicationIOError
Bases: OSError
OPC-UA communication I/O error.


exception OpcUaCommunicationTimeoutError
Bases: hvl_ccb.comm.opc.OpcUaCommunicationIOError
OPC-UA communication timeout error.

class OpcUaSubHandler
Bases: object
Base class for subscription handling of OPC events and data change events. Override methods from this class to add own handling capabilities.

To receive events from server for a subscription data_change and event methods are called directly from receiving thread. Do not do expensive, slow or network operation there. Create another thread if you need to do such a thing.

    datachange_notification(node, val, data)
    event_notification(event)
```

## hvl\_ccb.comm.serial

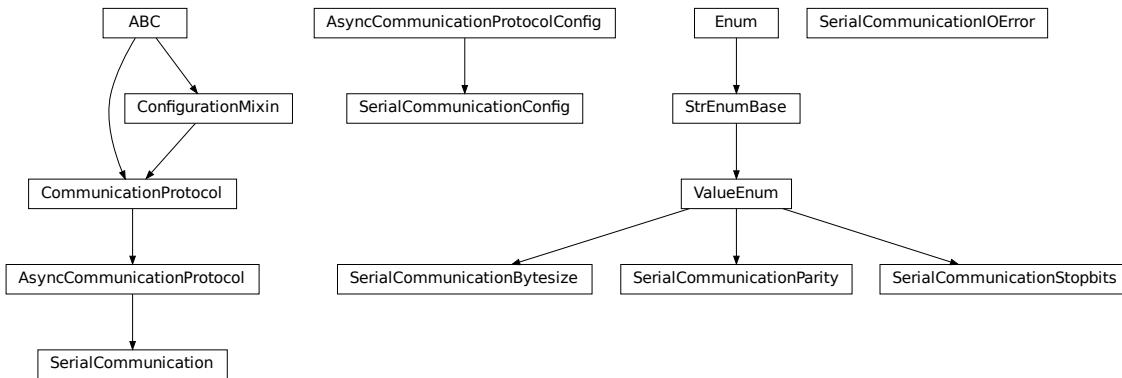

Communication protocol for serial ports. Makes use of the `pySerial` library.

**class** `SerialCommunication`(*configuration*)  
Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocol`

Implements the Communication Protocol for serial ports.

**close()**

Close the serial connection.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property is\_open**

Flag indicating if the serial port is open.

**Returns** `True` if the serial port is open, otherwise `False`

**open()**

Open the serial connection.

**Raises** `SerialCommunicationIOError` – when communication port cannot be opened.

**read\_bytes()** → bytes

Read the bytes from the serial port till the terminator is found. The input buffer may hold additional lines afterwards.

This method uses `self.access_lock` to ensure thread-safety.

**Returns** Bytes read from the serial port; `b''` if there was nothing to read.

**Raises** `SerialCommunicationIOError` – when communication port is not opened

**read\_single\_bytes(size: int = 1)** → bytes

Read the specified number of bytes from the serial port. The input buffer may hold additional data afterwards.

**Returns** Bytes read from the serial port; `b''` if there was nothing to read.

**write\_bytes(data: bytes)** → int

Write bytes to the serial port.

This method uses `self.access_lock` to ensure thread-safety.

**Parameters** `data` – data to write to the serial port

**Returns** number of bytes written

**Raises** `SerialCommunicationIOError` – when communication port is not opened

```
class SerialCommunicationBytesize(value=<object object>, names=None, module=None,
                                   type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Serial communication bytesize.

`EIGHTBITS` = 8

`FIVEBITS` = 5

`SEVENBITS` = 7

`SIXBITS` = 6

```
class SerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str =
                                'utf-8', encoding_error_handling: str = 'replace',
                                wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                default_n_attempts_read_text_nonempty: int = 10, port:
                                Optional[str] = None, baudrate: int = 9600, parity:
                                Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                Union[int, float, hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                Union[int, float] = 2)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

Configuration dataclass for `SerialCommunication`.

#### `Bytesize`

alias of `hvl_ccb.comm.serial.SerialCommunicationBytesize`

#### `Parity`

alias of `hvl_ccb.comm.serial.SerialCommunicationParity`

#### `Stopbits`

alias of `hvl_ccb.comm.serial.SerialCommunicationStopbits`

#### `baudrate: int = 9600`

Baudrate of the serial port

#### `bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8`

Size of a byte, 5 to 8

#### `clean_values()`

#### `create_serial_port() → serial.serialposix.Serial`

Create a serial port instance according to specification in this configuration

**Returns** Closed serial port instance

#### `force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.**parity:** Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'

Parity to be used for the connection.

**port:** Optional[str] = NonePort is a string referring to a COM-port (e.g. 'COM3') or a URL. The full list of capabilities is found [on the pyserial documentation](#).**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.**stopbits:** Union[int, float, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1

Stopbits setting, can be 1, 1.5 or 2.

**terminator\_str()** → str**timeout:** Union[int, float] = 2

Timeout in seconds for the serial port

**exception SerialCommunicationIOError**

Bases: OSError

Serial communication related I/O errors.

**class SerialCommunicationParity**(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)

Bases: hvl\_ccb.utils.enum.ValueEnum

Serial communication parity.

EVEN = 'E'

MARK = 'M'

NAMEs = { 'E': 'Even', 'M': 'Mark', 'N': 'None', 'O': 'Odd', 'S': 'Space' }

NONE = 'N'

ODD = 'O'

SPACE = 'S'


**class SerialCommunicationStopbits**(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)

Bases: hvl\_ccb.utils.enum.ValueEnum

Serial communication stopbits.

```
ONE = 1
ONE_POINT_FIVE = 1.5
TWO = 2
```

## hvl\_ccb.comm.telnet

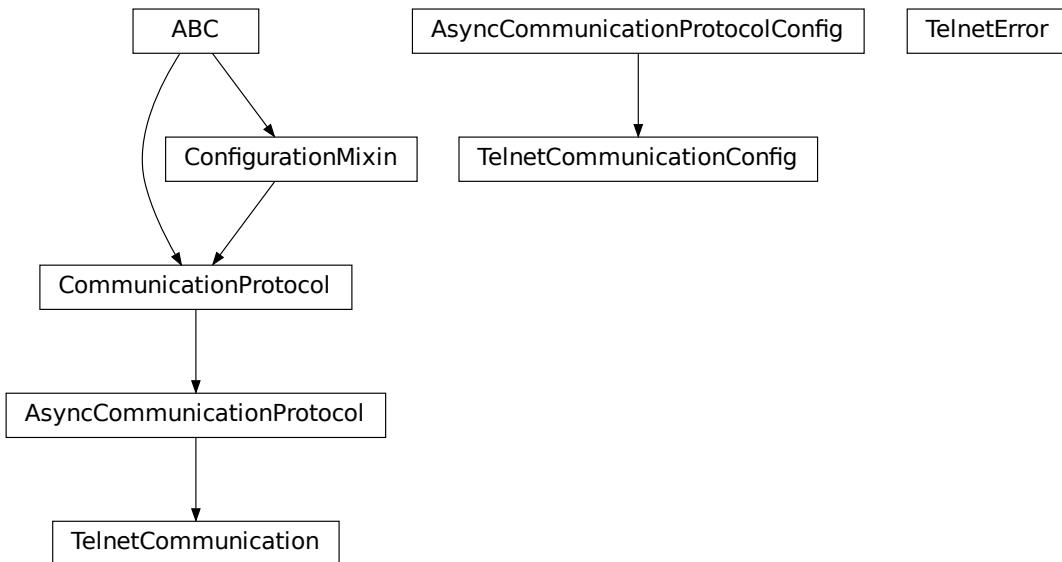

Communication protocol for telnet. Makes use of the [telnetlib](#) library.

```
class TelnetCommunication(configuration)
Bases: hvl_ccb.comm.base.AsyncCommunicationProtocol
```

Implements the Communication Protocol for telnet.

**close()**  
Close the telnet connection unless it is not closed.

**static config\_cls()**  
Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**property is\_open**  
Is the connection open?

**Returns** True for an open connection

**open()**  
Open the telnet connection unless it is not yet opened.

**read\_bytes() → bytes**  
Read data as *bytes* from the telnet connection.

**Returns** data from telnet connection

**Raises** `TelnetError` – when connection is not open, raises an Error during the communication

**write\_bytes** (`data: bytes`)

Write the data as `bytes` to the telnet connection.

**Parameters** `data` – Data to be sent.

**Raises** `TelnetError` – when connection is not open, raises an Error during the communication

```
class TelnetCommunicationConfig(terminator: bytes = b'\r\n', encoding: str =
    'utf-8', encoding_error_handling: str = 'replace',
    wait_sec_read_text_nonempty: Union[int, float] = 0.5,
    default_n_attempts_read_text_nonempty: int = 10, host:
    Optional[str] = None, port: int = 0, timeout: Union[int, float]
    = 0.2)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

Configuration dataclass for `TelnetCommunication`.

**clean\_values** ()

**create\_telnet** () → `Optional[telnetlib.Telnet]`

Create a telnet client :return: Opened Telnet object or None if connection is not possible

**force\_value** (`fieldname, value`)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**host:** `Optional[str] = None`

Host to connect to can be `localhost` or

**classmethod keys** () → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults** () → `Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port:** `int = 0`

Port at which the host is listening

**classmethod required\_keys** () → `Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**timeout:** `Union[int, float] = 0.2`


Timeout for reading a line

**exception TelnetError**

Bases: Exception

Telnet communication related errors.

**hvl\_ccb.comm.visa**

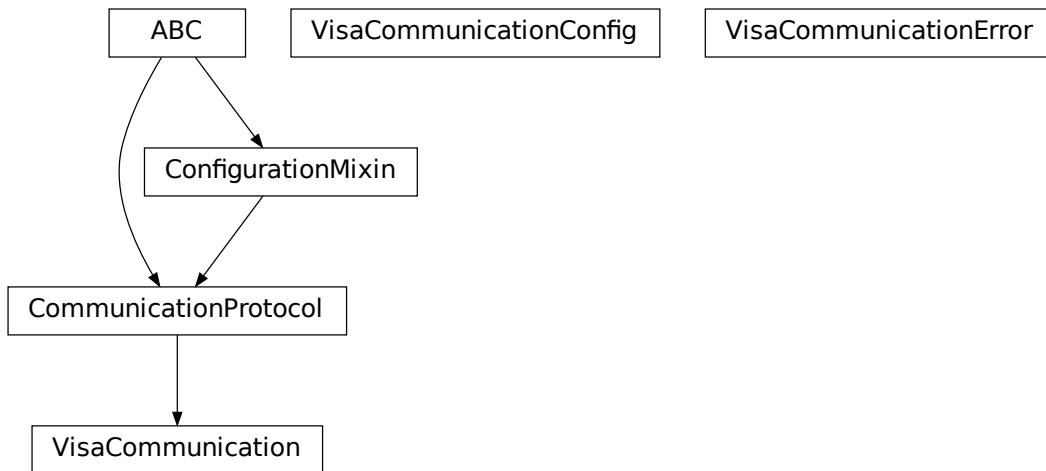

Communication protocol for VISA. Makes use of the pyvisa library. The backend can be NI-Visa or pyvisa-py.

Information on how to install a VISA backend can be found here: [https://pyvisa.readthedocs.io/en/master/getting\\_nivisa.html](https://pyvisa.readthedocs.io/en/master/getting_nivisa.html)

So far only TCPIP SOCKET and TCPIP INSTR interfaces are supported.

**class VisaCommunication(configuration)**

Bases: *hvl\_ccb.comm.base.CommunicationProtocol*

Implements the Communication Protocol for VISA / SCPI.

**MULTI\_COMMANDS\_MAX = 5**

The maximum of commands that can be sent in one round is 5 according to the VISA standard.

**MULTI\_COMMANDS\_SEPARATOR = ';'**

The character to separate two commands is ; according to the VISA standard.

**WAIT\_AFTER\_WRITE = 0.08**

Small pause in seconds to wait after write operations, allowing devices to really do what we tell them before continuing with further tasks.

**close() → None**

Close the VISA connection and invalidates the handle.

**static config\_cls() → Type[*hvl\_ccb.comm.visa.VisaCommunicationConfig*]**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**open () → None**  
 Open the VISA connection and create the resource.

**query (\*commands: str) → Union[str, Tuple[str, ...]]**  
 A combination of write(message) and read.

**Parameters** `commands` – list of commands

**Returns** list of values

**Raises** `VisaCommunicationError` – when connection was not started, or when trying to issue too many commands at once.

**spoll () → int**  
 Execute serial poll on the device. Reads the status byte register STB. This is a fast function that can be executed periodically in a polling fashion.

**Returns** integer representation of the status byte

**Raises** `VisaCommunicationError` – when connection was not started

**write (\*commands: str) → None**  
 Write commands. No answer is read or expected.

**Parameters** `commands` – one or more commands to send

**Raises** `VisaCommunicationError` – when connection was not started

---

**class VisaCommunicationConfig (host: str, interface\_type: Union[str, hvl\_ccb.comm.visa.VisaCommunicationConfig.InterfaceType], board: int = 0, port: int = 5025, timeout: int = 5000, chunk\_size: int = 204800, open\_timeout: int = 1000, write\_termination: str = '\n', read\_termination: str = '\n', visa\_backend: str = '')**

Bases: `object`

*VisaCommunication configuration dataclass.*

**class InterfaceType (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)**

Bases: `hvl_ccb.utils.enum.AutoNumberNameEnum`

Supported VISA Interface types.

**TCP/IP\_INSTR = 2**  
 VXI-11 protocol

**TCP/IP\_SOCKET = 1**  
 VISA-Raw protocol

**address (host: str, port: Optional[int] = None, board: Optional[int] = None) → str**  
 Address string specific to the VISA interface type.

**Parameters**

- `host` – host IP address
- `port` – optional TCP port
- `board` – optional board number

**Returns** address string

**property address**  
 Address string depending on the VISA protocol's configuration.

**Returns** address string corresponding to current configuration

**board: int = 0**  
 Board number is typically 0 and comes from old bus systems.

**chunk\_size: int = 204800**

Chunk size is the allocated memory for read operations. The standard is 20kB, and is increased per default here to 200kB. It is specified in bytes.

**clean\_values()****force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**host: str**

IP address of the VISA device. DNS names are currently unsupported.

**interface\_type: Union[str, hvl\_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]**  
Interface type of the VISA connection, being one of *InterfaceType*.**is\_configdataclass = True****classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**open\_timeout: int = 1000**

Timeout for opening the connection, in milli seconds.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 5025**

TCP port, standard is 5025.

**read\_termination: str = '\n'**

Read termination character.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**timeout: int = 5000**

Timeout for commands in milli seconds.

**visa\_backend: str = ''**

Specifies the path to the library to be used with PyVISA as a backend. Defaults to None, which is NI-VISA (if installed), or pyvisa-py (if NI-VISA is not found). To force the use of pyvisa-py, specify '@py' here.

**write\_termination: str = '\n'**

Write termination character.

**exception VisaCommunicationError**

Bases: Exception

Base class for VisaCommunication errors.

## Module contents

Communication protocols subpackage.


### **hvl\_ccb.dev**

#### Subpackages

##### **hvl\_ccb.dev.supercube**

#### Submodules

##### **hvl\_ccb.dev.supercube.base**

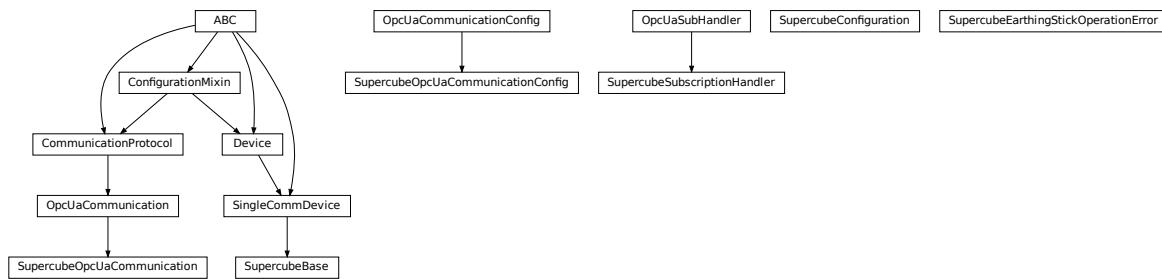

Base classes for the Supercube device.

**class SupercubeBase (com, dev\_config=None)**  
Bases: [hvl\\_ccb.dev.base.SingleCommDevice](#)

Base class for Supercube variants.

**static config\_cls()**  
Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**  
Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**display\_message\_board()** → None  
Display 15 newest messages

**display\_status\_board()** → None  
Display status board.

**get\_cee16\_socket()** → bool  
Read the on-state of the IEC CEE16 three-phase power socket.

**Returns** the on-state of the CEE16 power socket

**get\_door\_status** (*door: int*) → *hvl\_ccb.dev.supercube.constants.DoorStatus*  
Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

**Parameters** `door` – the door number (1..3)

**Returns** the door status

**get\_earthing\_rod\_status** (*earthing\_rod: int*) → *hvl\_ccb.dev.supercube.constants.EarthingRodStatus*  
Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

**Parameters** `earthing_rod` – the earthing rod number (1..3)

**Returns** the earthing rod status

**get\_earthing\_stick\_manual** (*number: int*) → *hvl\_ccb.dev.supercube.constants.EarthingStickOperation*  
Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** operation of the earthing stick in a manual operating mode (open == 0, close == 1)

**Raises** `ValueError` – when earthing stick number is not valid

**get\_earthing\_stick\_operating\_status** (*number: int*) → *hvl\_ccb.dev.supercube.constants.EarthingStickOperatingStatus*  
Get the operating status of an earthing stick.

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** earthing stick operating status (auto == 0, manual == 1)

**Raises** `ValueError` – when earthing stick number is not valid

**get\_earthing\_stick\_status** (*number: int*) → *hvl\_ccb.dev.supercube.constants.EarthingStickStatus*  
Get the status of an earthing stick, whether it is closed, open or undefined (moving).

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** earthing stick status

**Raises** `ValueError` – when earthing stick number is not valid

**get\_measurement\_ratio** (*channel: int*) → float  
Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

**Parameters** `channel` – number of the input channel (1..4)

**Returns** the ratio

**Raises** `ValueError` – when channel is not valid

**get\_measurement\_voltage** (*channel: int*) → float  
Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

**Parameters** `channel` – number of the input channel (1..4)

**Returns** measured voltage

**Raises** `ValueError` – when channel is not valid

**get\_status** () → *hvl\_ccb.dev.supercube.constants.SafetyStatus*  
Get the safety circuit status of the Supercube. :return: the safety status of the supercube's state machine.

**get\_support\_input** (*port: int, contact: int*) → bool

Get the state of a support socket input.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

**Returns** digital input read state

**Raises ValueError** – when port or contact number is not valid

**get\_support\_output** (*port: int, contact: int*) → bool

Get the state of a support socket output.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

**Returns** digital output read state

**Raises ValueError** – when port or contact number is not valid

**get\_t13\_socket** (*port: int*) → bool

Read the state of a SEV T13 power socket.

**Parameters** **port** – is the socket number, one of *constants.T13\_SOCKET\_PORTS*

**Returns** on-state of the power socket

**Raises ValueError** – when port is not valid

**operate** (*state: bool*) → None

Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

**Parameters** **state** – set operate state

**operate\_earthing\_stick** (*number: int, operation: hvl\_ccb.dev.supercube.constants.EarthingStickOperation*)

→ None

Operation of an earthing stick, which is set to manual operation. If an earthing stick is set to manual, it stays closed even if the system is in states RedReady or RedOperate.

**Parameters**

- **number** – number of the earthing stick (1..6)
- **operation** – earthing stick manual status (close or open)

**Raises SupercubeEarthingStickOperationError** – when operating status of given number's earthing stick is not manual

**quit\_error** () → None

Quits errors that are active on the Supercube.

**read** (*node\_id: str*)

Local wrapper for the OPC UA communication protocol read method.

**Parameters** **node\_id** – the id of the node to read.

**Returns** the value of the variable

**ready** (*state: bool*) → None

Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

**Parameters** **state** – set ready state

**set\_cee16\_socket** (*state: bool*) → None

Switch the IEC CEE16 three-phase power socket on or off.

**Parameters** **state** – desired on-state of the power socket

**Raises ValueError** – if state is not of type bool

**set\_message\_board** (*msgs: List[str], display\_board: bool = True*) → None

Fills messages into message board that display that 15 newest messages with a timestamp.

**Parameters**

- **msgs** – list of strings

- **display\_board** – display 15 newest messages if *True* (default)

**Raises ValueError** – if there are too many messages or the positions indices are invalid.

**set\_remote\_control** (*state: bool*) → None

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

**Parameters** **state** – desired remote control state

**set\_status\_board** (*msgs: List[str], pos: Optional[List[int]] = None, clear\_board: bool = True, display\_board: bool = True*) → None

Sets and displays a status board. The messages and the position of the message can be defined.

**Parameters**

- **msgs** – list of strings

- **pos** – list of integers [0...14]

- **clear\_board** – clear unspecified lines if *True* (default), keep otherwise

- **display\_board** – display new status board if *True* (default)

**Raises ValueError** – if there are too many messages or the positions indices are invalid.

**set\_support\_output** (*port: int, contact: int, state: bool*) → None

Set the state of a support output socket.

**Parameters**

- **port** – is the socket number (1..6)

- **contact** – is the contact on the socket (1..2)

- **state** – is the desired state of the support output

**Raises ValueError** – when port or contact number is not valid

**set\_support\_output\_impulse** (*port: int, contact: int, duration: float = 0.2, pos\_pulse: bool = True*) → None

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the *pos\_pulse* argument.

This function is blocking.

**Parameters**

- **port** – is the socket number (1..6)

- **contact** – is the contact on the socket (1..2)

- **duration** – is the length of the impulse in seconds

- **pos\_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

**Raises ValueError** – when port or contact number is not valid

**set\_t13\_socket** (*port: int, state: bool*) → None  
Set the state of a SEV T13 power socket.

#### Parameters

- **port** – is the socket number, one of *constants.T13\_SOCKET\_PORTS*
- **state** – is the desired on-state of the socket

**Raises ValueError** – when port is not valid or state is not of type bool

**start()** → None  
Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

**stop()** → None  
Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

**write** (*node\_id, value*) → None  
Local wrapper for the OPC UA communication protocol write method.

#### Parameters

- **node\_id** – the id of the node to read
- **value** – the value to write to the variable

**class SupercubeConfiguration** (*namespace\_index: int = 3, polling\_delay\_sec: Union[int, float] = 5.0, polling\_interval\_sec: Union[int, float] = 1.0*)

Bases: *object*

Configuration dataclass for the Supercube devices.

**clean\_values()**

**force\_value** (*fieldname, value*)  
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**namespace\_index: int = 3**  
Namespace of the OPC variables, typically this is 3 (coming from Siemens)

**classmethod optional\_defaults()** → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_delay\_sec: Union[int, float] = 5.0**

**polling\_interval\_sec: Union[int, float] = 1.0**

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception SupercubeEarthingStickOperationError**

Bases: Exception

**class SupercubeOpcUaCommunication(config)**

Bases: [hvl\\_ccb.comm.opc.OpcUaCommunication](#)

Communication protocol specification for Supercube devices.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class SupercubeOpcUaCommunicationConfig(host: str, endpoint\_name: str,**

**port: int = 4840, sub\_handler: hvl\_ccb.comm.opc.OpcUaSubHandler =**

**<hvl\_ccb.dev.supercube.base.SupercubeSubscriptionHandler**

**object>, update\_period: int = 500,**

**wait\_timeout\_retry\_sec: Union[int, float] = 1, max\_timeout\_retry\_nr: int = 5**

Bases: [hvl\\_ccb.comm.opc.OpcUaCommunicationConfig](#)

Communication protocol configuration for OPC UA, specifications for the Supercube devices.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field

- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**sub\_handler: hvl\_ccb.comm.opc.OpcUaSubHandler = <hvl\_ccb.dev.supercube.base.Supercube**

Subscription handler for data change events

**class SupercubeSubscriptionHandler**

Bases: [hvl\\_ccb.comm.opc.OpcUaSubHandler](#)

OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.


#### **datachange\_notification** (*node: opcua.common.node.Node, val, data*)

In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.

#### Parameters

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

### **hvl\_ccb.dev.supercube.constants**

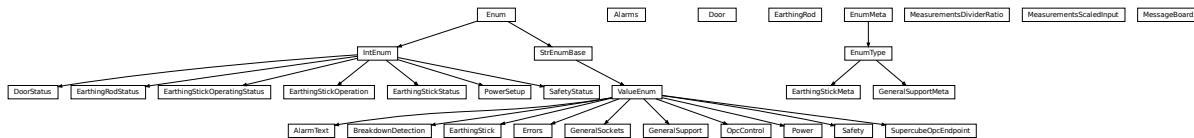

Constants, variable names for the Supercube OPC-connected devices.

```
class AlarmText (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the [AlarmText.get\(\)](#) method to retrieve the enum of an alarm number.

```

Alarm1 = 'STOP Emergency Stop 1'
Alarm10 = 'STOP Earthing stick 2 error while opening'
Alarm11 = 'STOP Earthing stick 3 error while opening'
Alarm12 = 'STOP Earthing stick 4 error while opening'
Alarm13 = 'STOP Earthing stick 5 error while opening'
Alarm14 = 'STOP Earthing stick 6 error while opening'
Alarm15 = 'STOP Earthing stick 1 error while closing'
Alarm16 = 'STOP Earthing stick 2 error while closing'
Alarm17 = 'STOP Earthing stick 3 error while closing'
Alarm18 = 'STOP Earthing stick 4 error while closing'
Alarm19 = 'STOP Earthing stick 5 error while closing'
Alarm2 = 'STOP Emergency Stop 2'
Alarm20 = 'STOP Earthing stick 6 error while closing'
Alarm21 = 'STOP Safety fence 1'
Alarm22 = 'STOP Safety fence 2'
Alarm23 = 'STOP OPC connection error'
Alarm24 = 'STOP Grid power failure'
```

```
Alarm25 = 'STOP UPS failure'
Alarm26 = 'STOP 24V PSU failure'
Alarm3 = 'STOP Emergency Stop 3'
Alarm4 = 'STOP Safety Switch 1 error'
Alarm41 = 'WARNING Door 1: Use earthing rod!'
Alarm42 = 'MESSAGE Door 1: Earthing rod is still in setup.'
Alarm43 = 'WARNING Door 2: Use earthing rod!'
Alarm44 = 'MESSAGE Door 2: Earthing rod is still in setup.'
Alarm45 = 'WARNING Door 3: Use earthing rod!'
Alarm46 = 'MESSAGE Door 3: Earthing rod is still in setup.'
Alarm47 = 'MESSAGE UPS charge < 85%'
Alarm48 = 'MESSAGE UPS running on battery'
Alarm5 = 'STOP Safety Switch 2 error'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error while opening'

classmethod get(alarm: int)
```

Get the attribute of this enum for an alarm number.

**Parameters** **alarm** – the alarm number

**Returns** the enum for the desired alarm number

```
not_defined = 'NO ALARM TEXT DEFINED'
```

```
class Alarms(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: hvl\_ccb.dev.supercube.constants.\_AlarmEnumBase

Alarms enumeration containing all variable NodeID strings for the alarm array.

```
Alarm1 = '"DB_Alarm_HMI"."Alarm1"'
Alarm10 = '"DB_Alarm_HMI"."Alarm10"'
Alarm100 = '"DB_Alarm_HMI"."Alarm100"'
Alarm101 = '"DB_Alarm_HMI"."Alarm101"'
Alarm102 = '"DB_Alarm_HMI"."Alarm102"'
Alarm103 = '"DB_Alarm_HMI"."Alarm103"'
Alarm104 = '"DB_Alarm_HMI"."Alarm104"'
Alarm105 = '"DB_Alarm_HMI"."Alarm105"'
Alarm106 = '"DB_Alarm_HMI"."Alarm106"'
Alarm107 = '"DB_Alarm_HMI"."Alarm107"'
Alarm108 = '"DB_Alarm_HMI"."Alarm108"'
```

```
Alarm109 = '"DB_Alarm_HMI"."Alarm109"'
Alarm11 = '"DB_Alarm_HMI"."Alarm11"'
Alarm110 = '"DB_Alarm_HMI"."Alarm110"'
Alarm111 = '"DB_Alarm_HMI"."Alarm111"'
Alarm112 = '"DB_Alarm_HMI"."Alarm112"'
Alarm113 = '"DB_Alarm_HMI"."Alarm113"'
Alarm114 = '"DB_Alarm_HMI"."Alarm114"'
Alarm115 = '"DB_Alarm_HMI"."Alarm115"'
Alarm116 = '"DB_Alarm_HMI"."Alarm116"'
Alarm117 = '"DB_Alarm_HMI"."Alarm117"'
Alarm118 = '"DB_Alarm_HMI"."Alarm118"'
Alarm119 = '"DB_Alarm_HMI"."Alarm119"'
Alarm12 = '"DB_Alarm_HMI"."Alarm12"'
Alarm120 = '"DB_Alarm_HMI"."Alarm120"'
Alarm121 = '"DB_Alarm_HMI"."Alarm121"'
Alarm122 = '"DB_Alarm_HMI"."Alarm122"'
Alarm123 = '"DB_Alarm_HMI"."Alarm123"'
Alarm124 = '"DB_Alarm_HMI"."Alarm124"'
Alarm125 = '"DB_Alarm_HMI"."Alarm125"'
Alarm126 = '"DB_Alarm_HMI"."Alarm126"'
Alarm127 = '"DB_Alarm_HMI"."Alarm127"'
Alarm128 = '"DB_Alarm_HMI"."Alarm128"'
Alarm129 = '"DB_Alarm_HMI"."Alarm129"'
Alarm13 = '"DB_Alarm_HMI"."Alarm13"'
Alarm130 = '"DB_Alarm_HMI"."Alarm130"'
Alarm131 = '"DB_Alarm_HMI"."Alarm131"'
Alarm132 = '"DB_Alarm_HMI"."Alarm132"'
Alarm133 = '"DB_Alarm_HMI"."Alarm133"'
Alarm134 = '"DB_Alarm_HMI"."Alarm134"'
Alarm135 = '"DB_Alarm_HMI"."Alarm135"'
Alarm136 = '"DB_Alarm_HMI"."Alarm136"'
Alarm137 = '"DB_Alarm_HMI"."Alarm137"'
Alarm138 = '"DB_Alarm_HMI"."Alarm138"'
Alarm139 = '"DB_Alarm_HMI"."Alarm139"'
Alarm14 = '"DB_Alarm_HMI"."Alarm14"'
Alarm140 = '"DB_Alarm_HMI"."Alarm140"'
```

```
Alarm141 = '"DB_Alarm_HMI"."Alarm141"'
Alarm142 = '"DB_Alarm_HMI"."Alarm142"'
Alarm143 = '"DB_Alarm_HMI"."Alarm143"'
Alarm144 = '"DB_Alarm_HMI"."Alarm144"'
Alarm145 = '"DB_Alarm_HMI"."Alarm145"'
Alarm146 = '"DB_Alarm_HMI"."Alarm146"'
Alarm147 = '"DB_Alarm_HMI"."Alarm147"'
Alarm148 = '"DB_Alarm_HMI"."Alarm148"'
Alarm149 = '"DB_Alarm_HMI"."Alarm149"'
Alarm15 = '"DB_Alarm_HMI"."Alarm15"'
Alarm150 = '"DB_Alarm_HMI"."Alarm150"'
Alarm151 = '"DB_Alarm_HMI"."Alarm151"'
Alarm16 = '"DB_Alarm_HMI"."Alarm16"'
Alarm17 = '"DB_Alarm_HMI"."Alarm17"'
Alarm18 = '"DB_Alarm_HMI"."Alarm18"'
Alarm19 = '"DB_Alarm_HMI"."Alarm19"'
Alarm2 = '"DB_Alarm_HMI"."Alarm2"'
Alarm20 = '"DB_Alarm_HMI"."Alarm20"'
Alarm21 = '"DB_Alarm_HMI"."Alarm21"'
Alarm22 = '"DB_Alarm_HMI"."Alarm22"'
Alarm23 = '"DB_Alarm_HMI"."Alarm23"'
Alarm24 = '"DB_Alarm_HMI"."Alarm24"'
Alarm25 = '"DB_Alarm_HMI"."Alarm25"'
Alarm26 = '"DB_Alarm_HMI"."Alarm26"'
Alarm27 = '"DB_Alarm_HMI"."Alarm27"'
Alarm28 = '"DB_Alarm_HMI"."Alarm28"'
Alarm29 = '"DB_Alarm_HMI"."Alarm29"'
Alarm3 = '"DB_Alarm_HMI"."Alarm3"'
Alarm30 = '"DB_Alarm_HMI"."Alarm30"'
Alarm31 = '"DB_Alarm_HMI"."Alarm31"'
Alarm32 = '"DB_Alarm_HMI"."Alarm32"'
Alarm33 = '"DB_Alarm_HMI"."Alarm33"'
Alarm34 = '"DB_Alarm_HMI"."Alarm34"'
Alarm35 = '"DB_Alarm_HMI"."Alarm35"'
Alarm36 = '"DB_Alarm_HMI"."Alarm36"'
Alarm37 = '"DB_Alarm_HMI"."Alarm37"'
```

```
Alarm38 = '"DB_Alarm_HMI"."Alarm38"'
Alarm39 = '"DB_Alarm_HMI"."Alarm39"'
Alarm4 = '"DB_Alarm_HMI"."Alarm4"'
Alarm40 = '"DB_Alarm_HMI"."Alarm40"'
Alarm41 = '"DB_Alarm_HMI"."Alarm41"'
Alarm42 = '"DB_Alarm_HMI"."Alarm42"'
Alarm43 = '"DB_Alarm_HMI"."Alarm43"'
Alarm44 = '"DB_Alarm_HMI"."Alarm44"'
Alarm45 = '"DB_Alarm_HMI"."Alarm45"'
Alarm46 = '"DB_Alarm_HMI"."Alarm46"'
Alarm47 = '"DB_Alarm_HMI"."Alarm47"'
Alarm48 = '"DB_Alarm_HMI"."Alarm48"'
Alarm49 = '"DB_Alarm_HMI"."Alarm49"'
Alarm5 = '"DB_Alarm_HMI"."Alarm5"'
Alarm50 = '"DB_Alarm_HMI"."Alarm50"'
Alarm51 = '"DB_Alarm_HMI"."Alarm51"'
Alarm52 = '"DB_Alarm_HMI"."Alarm52"'
Alarm53 = '"DB_Alarm_HMI"."Alarm53"'
Alarm54 = '"DB_Alarm_HMI"."Alarm54"'
Alarm55 = '"DB_Alarm_HMI"."Alarm55"'
Alarm56 = '"DB_Alarm_HMI"."Alarm56"'
Alarm57 = '"DB_Alarm_HMI"."Alarm57"'
Alarm58 = '"DB_Alarm_HMI"."Alarm58"'
Alarm59 = '"DB_Alarm_HMI"."Alarm59"'
Alarm6 = '"DB_Alarm_HMI"."Alarm6"'
Alarm60 = '"DB_Alarm_HMI"."Alarm60"'
Alarm61 = '"DB_Alarm_HMI"."Alarm61"'
Alarm62 = '"DB_Alarm_HMI"."Alarm62"'
Alarm63 = '"DB_Alarm_HMI"."Alarm63"'
Alarm64 = '"DB_Alarm_HMI"."Alarm64"'
Alarm65 = '"DB_Alarm_HMI"."Alarm65"'
Alarm66 = '"DB_Alarm_HMI"."Alarm66"'
Alarm67 = '"DB_Alarm_HMI"."Alarm67"'
Alarm68 = '"DB_Alarm_HMI"."Alarm68"'
Alarm69 = '"DB_Alarm_HMI"."Alarm69"'
Alarm7 = '"DB_Alarm_HMI"."Alarm7"'
```

```
Alarm70 = '"DB_Alarm_HMI"."Alarm70"'
Alarm71 = '"DB_Alarm_HMI"."Alarm71"'
Alarm72 = '"DB_Alarm_HMI"."Alarm72"'
Alarm73 = '"DB_Alarm_HMI"."Alarm73"'
Alarm74 = '"DB_Alarm_HMI"."Alarm74"'
Alarm75 = '"DB_Alarm_HMI"."Alarm75"'
Alarm76 = '"DB_Alarm_HMI"."Alarm76"'
Alarm77 = '"DB_Alarm_HMI"."Alarm77"'
Alarm78 = '"DB_Alarm_HMI"."Alarm78"'
Alarm79 = '"DB_Alarm_HMI"."Alarm79"'
Alarm8 = '"DB_Alarm_HMI"."Alarm8"'
Alarm80 = '"DB_Alarm_HMI"."Alarm80"'
Alarm81 = '"DB_Alarm_HMI"."Alarm81"'
Alarm82 = '"DB_Alarm_HMI"."Alarm82"'
Alarm83 = '"DB_Alarm_HMI"."Alarm83"'
Alarm84 = '"DB_Alarm_HMI"."Alarm84"'
Alarm85 = '"DB_Alarm_HMI"."Alarm85"'
Alarm86 = '"DB_Alarm_HMI"."Alarm86"'
Alarm87 = '"DB_Alarm_HMI"."Alarm87"'
Alarm88 = '"DB_Alarm_HMI"."Alarm88"'
Alarm89 = '"DB_Alarm_HMI"."Alarm89"'
Alarm9 = '"DB_Alarm_HMI"."Alarm9"'
Alarm90 = '"DB_Alarm_HMI"."Alarm90"'
Alarm91 = '"DB_Alarm_HMI"."Alarm91"'
Alarm92 = '"DB_Alarm_HMI"."Alarm92"'
Alarm93 = '"DB_Alarm_HMI"."Alarm93"'
Alarm94 = '"DB_Alarm_HMI"."Alarm94"'
Alarm95 = '"DB_Alarm_HMI"."Alarm95"'
Alarm96 = '"DB_Alarm_HMI"."Alarm96"'
Alarm97 = '"DB_Alarm_HMI"."Alarm97"'
Alarm98 = '"DB_Alarm_HMI"."Alarm98"'
Alarm99 = '"DB_Alarm_HMI"."Alarm99"'

class BreakdownDetection(value=<object object>, names=None, module=None, type=None,
                        start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Node ID strings for the breakdown detection.

TODO: these variable NodeIDs are not tested and/or correct yet.

```

activated = '"Ix_Allg_Breakdown_activated"'
Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the system or not.

reset = '"Qx_Allg_Breakdown_reset"'
Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

triggered = '"Ix_Allg_Breakdown_triggered"'
Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using the safety circuit state, therefore no method is implemented to read this out directly.

class Door(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._DoorEnumBase

Variable NodeID strings for doors.

status_1 = '"DB_Safety_Circuit"."Door_1"."si_HMI_status"'
status_2 = '"DB_Safety_Circuit"."Door_2"."si_HMI_status"'
status_3 = '"DB_Safety_Circuit"."Door_3"."si_HMI_status"'

class DoorStatus(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for doors.

closed = 2
Door is closed, but not locked.

error = 4
Door has an error or was opened in locked state (either with emergency stop or from the inside).

inactive = 0
not enabled in Supercube HMI setup, this door is not supervised.

locked = 3
Door is closed and locked (safe state).

open = 1
Door is open.

class EarthingRod(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._DoorEnumBase

Variable NodeID strings for earthing rods.

status_1 = '"DB_Safety_Circuit"."Door_1"."Ix_earthingrod"'
status_2 = '"DB_Safety_Circuit"."Door_2"."Ix_earthingrod"'
status_3 = '"DB_Safety_Circuit"."Door_3"."Ix_earthingrod"'

class EarthingRodStatus(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for earthing rods.

experiment_blocked = 0
earthing rod is somewhere in the experiment and blocks the start of the experiment

experiment_ready = 1
earthing rod is hanging next to the door, experiment is ready to operate

```

```
class EarthingStick(value=<object object>, names=None, module=None, type=None, start=1,
                      boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum
```

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the earthing in manual mode.

```
classmethod manual(number: int)
```

Get the manual enum instance for an earthing stick number.

**Parameters** **number** – the earthing stick (1..6)

**Returns** the manual instance

**Raises** **ValueError** – when earthing stick number is not valid

```
manual_1 = '"DB_Safety_Circuit"."Earthstick_1"."sx_earthing_manually"'
manual_2 = '"DB_Safety_Circuit"."Earthstick_2"."sx_earthing_manually"'
manual_3 = '"DB_Safety_Circuit"."Earthstick_3"."sx_earthing_manually"'
manual_4 = '"DB_Safety_Circuit"."Earthstick_4"."sx_earthing_manually"'
manual_5 = '"DB_Safety_Circuit"."Earthstick_5"."sx_earthing_manually"'
manual_6 = '"DB_Safety_Circuit"."Earthstick_6"."sx_earthing_manually"'
classmethod manuals() → Tuple[hvl\_ccb.dev.supercube.constants.EarthingStick, ...]
```

Get all earthing stick manual instances.

**Returns** tuple of manual instances

```
property number
```

Get corresponding earthing stick number.

**Returns** earthing stick number (1..6)

```
classmethod operating_status(number: int)
```

Get the operating status enum instance for an earthing stick number.

**Parameters** **number** – the earthing stick (1..6)

**Returns** the operating status instance

**Raises** **ValueError** – when earthing stick number is not valid

```
operating_status_1 = '"DB_Safety_Circuit"."Earthstick_1"."sx_manual_control_active"'
operating_status_2 = '"DB_Safety_Circuit"."Earthstick_2"."sx_manual_control_active"'
operating_status_3 = '"DB_Safety_Circuit"."Earthstick_3"."sx_manual_control_active"'
operating_status_4 = '"DB_Safety_Circuit"."Earthstick_4"."sx_manual_control_active"'
operating_status_5 = '"DB_Safety_Circuit"."Earthstick_5"."sx_manual_control_active"'
operating_status_6 = '"DB_Safety_Circuit"."Earthstick_6"."sx_manual_control_active"'
classmethod operating_statuses() → Tuple[hvl\_ccb.dev.supercube.constants.EarthingStick,
                                              ...]
```

Get all earthing stick operating status instances.

**Returns** tuple of operating status instances

```
classmethod range() → Sequence[int]
```

Integer range of all earthing sticks.

**Returns** sequence of earthing sticks numbers

---

**classmethod** **status** (*number: int*) → *hvl\_ccb.dev.supercube.constants.EarthingStick*  
Get the status enum instance for an earthing stick number.

**Parameters** **number** – the earthing stick (1..6)

**Returns** the status instance

**Raises** **ValueError** – when earthing stick number is not valid

```
status_1 = '"DB_Safety_Circuit"."Earthstick_1"."si_HMI_Status"'
status_2 = '"DB_Safety_Circuit"."Earthstick_2"."si_HMI_Status"'
status_3 = '"DB_Safety_Circuit"."Earthstick_3"."si_HMI_Status"'
status_4 = '"DB_Safety_Circuit"."Earthstick_4"."si_HMI_Status"'
status_5 = '"DB_Safety_Circuit"."Earthstick_5"."si_HMI_Status"'
status_6 = '"DB_Safety_Circuit"."Earthstick_6"."si_HMI_Status"'
```

**classmethod** **statuses** () → Tuple[*hvl\_ccb.dev.supercube.constants.EarthingStick*, ...]

Get all earthing stick status instances.

**Returns** tuple of status instances

**class** **EarthingStickMeta** (*clsname, bases, clsdict, \*\*kwargs*)

Bases: aenum.EnumType

**class** **EarthingStickOperatingStatus** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: aenum.IntEnum

Operating Status for an earthing stick. Stick can be used in auto or manual mode.

**auto** = 0

**manual** = 1

**class** **EarthingStickOperation** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: aenum.IntEnum

Operation of the earthing stick in manual operating mode. Can be closed or opened.

**close** = 1

**open** = 0

**class** **EarthingStickStatus** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: aenum.IntEnum

Status of an earthing stick. These are the possible values in the status integer e.g. in *EarthingStick.status\_1*.

**closed** = 1

Earthing is closed (safe).

**error** = 3

Earthing is in error, e.g. when the stick did not close correctly or could not open.

**inactive** = 0

Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be enabled in the Supercube HMI setup.

```
open = 2
    Earthing is open (not safe).

class Errors(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Variable NodeID strings for information regarding error, warning and message handling.

message = '"DB_Message_Buffer"."Info_active"'
    Boolean read-only variable telling if a message is active.

quit = '"DB_Message_Buffer"."Reset_button"'
    Writable boolean for the error quit button.

stop = '"DB_Message_Buffer"."Stop_active"'
    Boolean read-only variable telling if a stop is active.

warning = '"DB_Message_Buffer"."Warning_active"'
    Boolean read-only variable telling if a warning is active.

class GeneralSockets(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

NodeID strings for the power sockets (3x T13 and 1xCEE16).

cee16 = '"Qx_Allg_Socket_CEE16"'
    CEE16 socket (writeable boolean).

t13_1 = '"Qx_Allg_Socket_T13_1"'
    SEV T13 socket No. 1 (writable boolean).

t13_2 = '"Qx_Allg_Socket_T13_2"'
    SEV T13 socket No. 2 (writable boolean).

t13_3 = '"Qx_Allg_Socket_T13_3"'
    SEV T13 socket No. 3 (writable boolean).

class GeneralSupport(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

NodeID strings for the support inputs and outputs.

classmethod contact_range() → Sequence[int]
    Integer range of all contacts.

        Returns sequence of contact numbers

in_1_1 = '"Ix_Allg_Support1_1"'
in_1_2 = '"Ix_Allg_Support1_2"'
in_2_1 = '"Ix_Allg_Support2_1"'
in_2_2 = '"Ix_Allg_Support2_2"'
in_3_1 = '"Ix_Allg_Support3_1"'
in_3_2 = '"Ix_Allg_Support3_2"'
in_4_1 = '"Ix_Allg_Support4_1"'
in_4_2 = '"Ix_Allg_Support4_2"'
in_5_1 = '"Ix_Allg_Support5_1"'
```

```

in_5_2 = '"Ix_Allg_Support5_2"'
in_6_1 = '"Ix_Allg_Support6_1"'
in_6_2 = '"Ix_Allg_Support6_2"'

```

**classmethod input** (*port: int, contact: int*)  
Get the NodeID string for a support input.

#### Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

**Returns** the node id string

**Raises ValueError** – when port or contact number is not valid

```

out_1_1 = "Qx_Allg_Support1_1"
out_1_2 = "Qx_Allg_Support1_2"
out_2_1 = "Qx_Allg_Support2_1"
out_2_2 = "Qx_Allg_Support2_2"
out_3_1 = "Qx_Allg_Support3_1"
out_3_2 = "Qx_Allg_Support3_2"
out_4_1 = "Qx_Allg_Support4_1"
out_4_2 = "Qx_Allg_Support4_2"
out_5_1 = "Qx_Allg_Support5_1"
out_5_2 = "Qx_Allg_Support5_2"
out_6_1 = "Qx_Allg_Support6_1"
out_6_2 = "Qx_Allg_Support6_2"

```

**classmethod output** (*port: int, contact: int*)  
Get the NodeID string for a support output.

#### Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

**Returns** the node id string

**Raises ValueError** – when port or contact number is not valid

**classmethod port\_range()** → Sequence[int]

Integer range of all ports.

**Returns** sequence of port numbers

**class GeneralSupportMeta** (*clsname, bases, clsdict, \*\*kwargs*)

Bases: aenum.EnumType

**class MeasurementsDividerRatio** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: hvl\_ccb.dev.supercube.constants.\_InputEnumBase

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

```
input_1 = '"DB_Measurements"."si_Divider_Ratio_1"'
input_2 = '"DB_Measurements"."si_Divider_Ratio_2"'
input_3 = '"DB_Measurements"."si_Divider_Ratio_3"'
input_4 = '"DB_Measurements"."si_Divider_Ratio_4"'

class MeasurementsScaledInput (value=<object object>, names=None, module=None, type=None,
                               start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._InputEnumBase
```

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in *MeasurementsDividerRatio*.

```
input_1 = '"DB_Measurements"."si_scaled_Voltage_Input_1"'
input_2 = '"DB_Measurements"."si_scaled_Voltage_Input_2"'
input_3 = '"DB_Measurements"."si_scaled_Voltage_Input_3"'
input_4 = '"DB_Measurements"."si_scaled_Voltage_Input_4"'

class MessageBoard (value=<object object>, names=None, module=None, type=None, start=1, bound-
                     ary=None)
Bases: hvl_ccb.dev.supercube.constants._LineEnumBase
```

Variable NodeID strings for message board lines.

```
line_1 = '"DB_OP_C_Connection"."Is_status_Line_1"'
line_10 = '"DB_OP_C_Connection"."Is_status_Line_10"'
line_11 = '"DB_OP_C_Connection"."Is_status_Line_11"'
line_12 = '"DB_OP_C_Connection"."Is_status_Line_12"'
line_13 = '"DB_OP_C_Connection"."Is_status_Line_13"'
line_14 = '"DB_OP_C_Connection"."Is_status_Line_14"'
line_15 = '"DB_OP_C_Connection"."Is_status_Line_15"'
line_2 = '"DB_OP_C_Connection"."Is_status_Line_2"'
line_3 = '"DB_OP_C_Connection"."Is_status_Line_3"'
line_4 = '"DB_OP_C_Connection"."Is_status_Line_4"'
line_5 = '"DB_OP_C_Connection"."Is_status_Line_5"'
line_6 = '"DB_OP_C_Connection"."Is_status_Line_6"'
line_7 = '"DB_OP_C_Connection"."Is_status_Line_7"'
line_8 = '"DB_OP_C_Connection"."Is_status_Line_8"'
line_9 = '"DB_OP_C_Connection"."Is_status_Line_9"'
```

```
class OpcControl (value=<object object>, names=None, module=None, type=None, start=1, bound-
                      ary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Variable NodeID strings for supervision of the OPC connection from the controlling workstation to the Super-cube.

```
active = '"DB_OP_C_Connection"."sx_OP_C_active"'
writable boolean to enable OPC remote control and display a message window on the Supercube HMI.
```

```

live = '"DB_OPC_Connection".sx_OPC_lifebit"'

class Power(value=<object object>, names=None, module=None, type=None, start=1, bound-
ary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Variable NodeID strings concerning power data.

TODO: these variable NodeIDs are not tested and/or correct yet, they don't exist yet on Supercube side.

current_primary = 'Qr_Power_FU_actual_Current'
    Primary current in ampere, measured by the frequency converter. (read-only)

frequency = 'Ir_Power_FU_Frequency'
    Frequency converter output frequency. (read-only)

setup = 'Qi_Power_Setup'
    Power setup that is configured using the Supercube HMI. The value corresponds to the ones in
    PowerSetup. (read-only)

voltage_max = 'Iw_Power_max_Voltage'
    Maximum voltage allowed by the current experimental setup. (read-only)

voltage_primary = 'Qr_Power_FU_actual_Voltage'
    Primary voltage in volts, measured by the frequency converter at its output. (read-only)

voltage_slope = 'Ir_Power_dUdt'
    Voltage slope in V/s.

voltage_target = 'Ir_Power_Target_Voltage'
    Target voltage setpoint in V.

class PowerSetup(value=<object object>, names=None, module=None, type=None, start=1, bound-
ary=None)
Bases: aenum.IntEnum

Possible power setups corresponding to the value of variable Power.setup.

AC_DoubleStage_150kV = 4
    AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum
    voltage of 150kV.

AC_DoubleStage_200kV = 5
    AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

AC_SingleStage_100kV = 3
    AC voltage with MWB transformer set to 100kV maximum voltage.

AC_SingleStage_50kV = 2
    AC voltage with MWB transformer set to 50kV maximum voltage.

DC_DoubleStage_280kV = 8
    DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single
    stage transformer with Greinacher voltage doubling rectifier)

DC_SingleStage_140kV = 7
    DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

External = 1
    External power supply fed through blue CEE32 input using isolation transformer and safety switches of
    the Supercube, or using an external safety switch attached to the Supercube Type B.

```

**Internal = 6**

Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

**NoPower = 0**

No safety switches, use only safety components (doors, fence, earthing...) without any power.

**class Safety**(*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl\_ccb.utils.enum.ValueEnum*

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

**status = '"DB\_Safety\_Circuit"."si\_safe\_status"**

Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in *SafetyStatus*.

**switch\_to\_operate = '"DB\_Safety\_Circuit"."sx\_safe\_switch\_to\_operate"**

Writable boolean for switching to Red Operate (locket, HV on) state.

**switch\_to\_ready = '"DB\_Safety\_Circuit"."sx\_safe\_switch\_to\_ready"**

Writable boolean for switching to Red Ready (locked, HV off) state.

**class SafetyStatus**(*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *aenum.IntEnum*

Safety status values that are possible states returned from *hvl\_ccb.dev.supercube.base.Supercube.get\_status()*. These values correspond to the states of the Supercube’s safety circuit statemachine.

**Error = 6**

System is in error mode.

**GreenNotReady = 1**

System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

**GreenReady = 2**

System is safe and all safety elements are in place to be able to switch to *ready*.

**Initializing = 0**

System is initializing or booting.

**QuickStop = 5**

Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

**RedOperate = 4**

System is locked in red state and in *operate* mode, i.e. high voltage on.

**RedReady = 3**

System is locked in red state and *ready* to go to *operate* mode.

**class SupercubeOpcEndpoint**(*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)


Bases: *hvl\_ccb.utils.enum.ValueEnum*

OPC Server Endpoint strings for the supercube variants.

**A = 'Supercube Typ A'****B = 'Supercube Typ B'**

```
T13_SOCKET_PORTS = (1, 2, 3)
Port numbers of SEV T13 power socket
```

### `hvl_ccb.dev.supercube.typ_a`

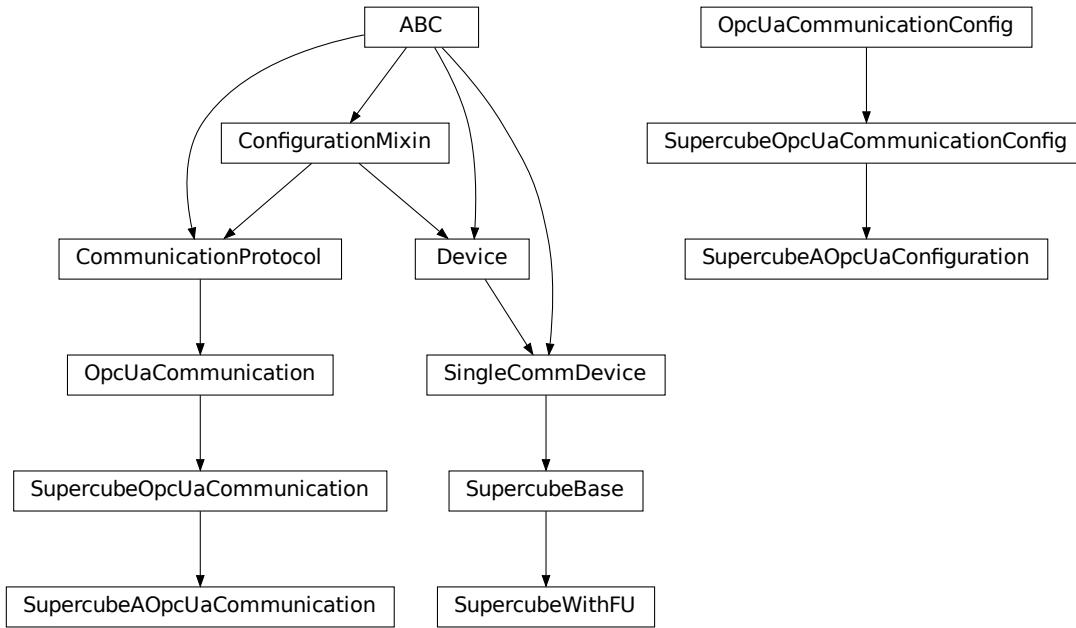

Supercube Typ A module.

```
class SupercubeAOpcaCommunication(config)
Bases: hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunication

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

class SupercubeAOpcaConfiguration(host: str, endpoint_name: str = 'Supercube Typ A', port: int = 4840, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object at 0x7f37219f9590>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)
Bases: hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunicationConfig

endpoint_name: str = 'Supercube Typ A'
    Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.
```

### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**class SupercubeWithFU**(com, dev\_config=None)

Bases: *hvl\_ccb.dev.supercube.base.SuperCubeBase*

Variant A of the Supercube with frequency converter.

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**fso\_reset()** → None

**TODO:** test fso\_reset with device

Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.

**get\_frequency()** → float

**TODO:** test get\_frequency with device

Read the electrical frequency of the current Supercube setup.

**Returns** the frequency in Hz

**get\_fso\_active()** → bool

**TODO:** test get\_fso\_active with device

Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.

**Returns** state of the FSO functionality

**get\_max\_voltage()** → float

**TODO:** test get\_max\_voltage with device

Reads the maximum voltage of the setup and returns in V.

**Returns** the maximum voltage of the setup in V.

**get\_power\_setup()** → *hvl\_ccb.dev.supercube.constants.PowerSetup*

**TODO:** test get\_power\_setup with device

Return the power setup selected in the Supercube's settings.

**Returns** the power setup

`get_primary_current () → float`

**TODO: get\_primary\_current with device**

Read the current primary current at the output of the frequency converter ( before transformer).

**Returns** primary current in A

`get_primary_voltage () → float`

**TODO: test get\_primary\_voltage with device**

Read the current primary voltage at the output of the frequency converter ( before transformer).

**Returns** primary voltage in V

`get_target_voltage () → float`

**TODO: test get\_target\_voltage with device**

Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to `set_target_voltage ()`.

**Returns** the setpoint voltage in V.

`set_slope (slope: float) → None`

**TODO: test set\_slope with device**

Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.

**Parameters** `slope` – voltage slope in V/s (0..15)


`set_target_voltage (volt_v: float) → None`

**TODO: test set\_target\_voltage with device**

Set the output voltage to a defined value in V.

**Parameters** `volt_v` – the desired voltage in V

## `hvl_ccb.dev.supercube.typ_b`

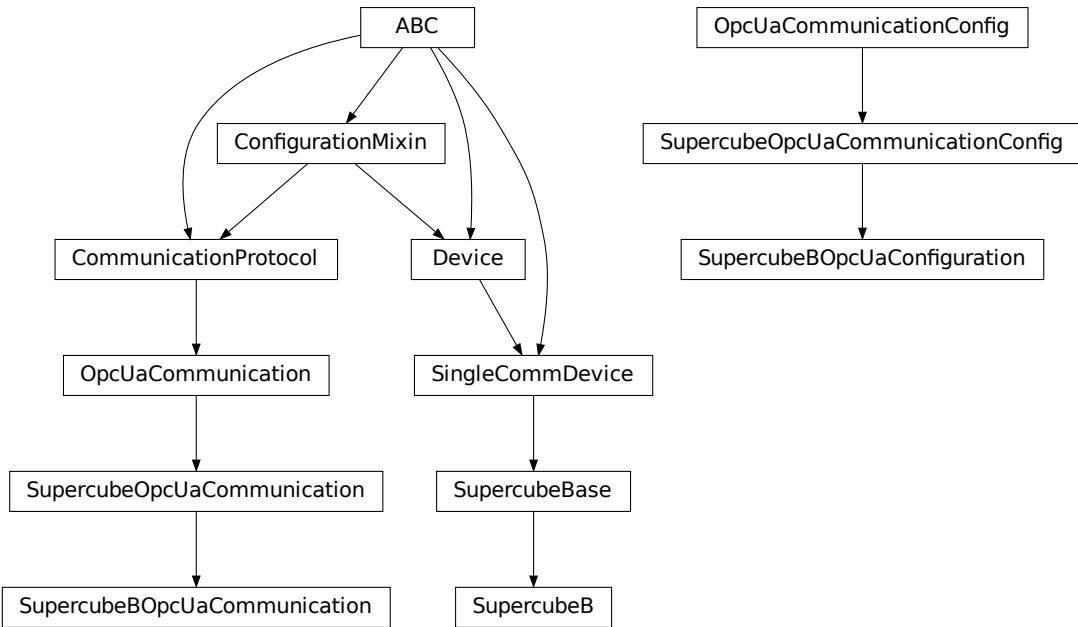

Supercube Typ B module.

```
class SupercubeB(com, dev_config=None)
Bases: hvl_ccb.dev.supercube.base.SupercubeBase
```

Variant B of the Supercube without frequency converter but external safety switches.

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

```
class SupercubeBOpcUaCommunication(config)
```

```
Bases: hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunication
```

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class SupercubeBOpcUaConfiguration(host: str, endpoint_name: str = 'Supercube Typ B', port: int = 4840, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object at 0x7f37219f9590>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)
```

```
Bases: hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunicationConfig
```

```
endpoint_name: str = 'Supercube Typ B'
```

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

**force\_value** (*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys () → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults () → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.


**classmethod required\_keys () → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**Module contents**

Supercube package with implementation for system versions from 2019 on (new concept with hard-PLC Siemens S7-1500 as CPU).

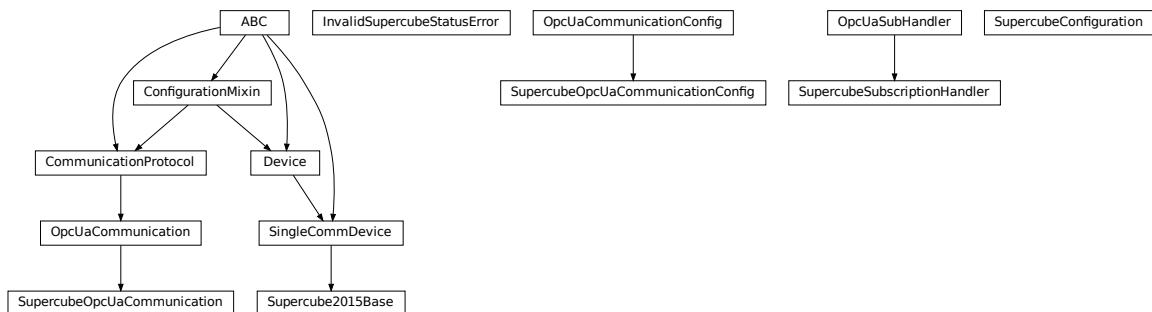**hvl\_ccb.dev.supercube2015****Submodules****hvl\_ccb.dev.supercube2015.base**

Base classes for the Supercube device.

**exception InvalidSupercubeStatusError**

Bases: Exception

Exception raised when supercube has invalid status.

**class Supercube2015Base (com, dev\_config=None)**

Bases: *hvl\_ccb.dev.base.SingleCommDevice*

Base class for Supercube variants.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**get\_cee16\_socket () → bool**

Read the on-state of the IEC CEE16 three-phase power socket.

**Returns** the on-state of the CEE16 power socket

**get\_door\_status (door: int) → hvl\_ccb.dev.supercube2015.constants.DoorStatus**

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

**Parameters** `door` – the door number (1..3)

**Returns** the door status

**get\_earthing\_manual (number: int) → bool**

Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** earthing stick manual status

**get\_earthing\_status (number: int) → int**

Get the status of an earthing stick, whether it is closed, open or undefined (moving).

**Parameters** `number` – number of the earthing stick (1..6)

**Returns** earthing stick status; see `constants.EarthingStickStatus`

**get\_measurement\_ratio (channel: int) → float**

Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

**Attention:** Supercube 2015 does not have a separate ratio for every analog input. Therefore there is only one ratio for `channel = 1`.

**Parameters** `channel` – number of the input channel (1..4)

**Returns** the ratio

**get\_measurement\_voltage (channel: int) → float**

Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

**Attention:** In contrast to the *new* Supercube, the old one returns here the input voltage read at the ADC. It is not scaled by a factor.

**Parameters** `channel` – number of the input channel (1..4)

**Returns** measured voltage

**get\_status ()** → int

Get the safety circuit status of the Supercube.

**Returns** the safety status of the supercube's state machine; see *constants.SafetyStatus*.

**get\_support\_input (port: int, contact: int)** → bool

Get the state of a support socket input.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

**Returns** digital input read state

**get\_support\_output (port: int, contact: int)** → bool

Get the state of a support socket output.

**Parameters**

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

**Returns** digital output read state

**get\_t13\_socket (port: int)** → bool

Read the state of a SEV T13 power socket.

**Parameters** **port** – is the socket number, one of *constants.T13\_SOCKET\_PORTS*

**Returns** on-state of the power socket

**horn (state: bool)** → None

Turns acoustic horn on or off.

**Parameters** **state** – Turns horn on (True) or off (False)

**operate (state: bool)** → None

Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

**Parameters** **state** – set operate state

**quit\_error ()** → None

Quits errors that are active on the Supercube.

**read (node\_id: str)**

Local wrapper for the OPC UA communication protocol read method.

**Parameters** **node\_id** – the id of the node to read.

**Returns** the value of the variable

**ready (state: bool)** → None

Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

**Parameters** **state** – set ready state

**set\_cee16\_socket (state: bool)** → None

Switch the IEC CEE16 three-phase power socket on or off.

**Parameters** **state** – desired on-state of the power socket

**Raises** **ValueError** – if state is not of type bool

**set\_earthing\_manual** (*number: int, manual: bool*) → None  
Set the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

### Parameters

- **number** – number of the earthing stick (1..6)
- **manual** – earthing stick manual status (True or False)

**set\_remote\_control** (*state: bool*) → None

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

### Parameters **state** – desired remote control state

**set\_support\_output** (*port: int, contact: int, state: bool*) → None

Set the state of a support output socket.

### Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **state** – is the desired state of the support output

**set\_support\_output\_impulse** (*port: int, contact: int, duration: float = 0.2, pos\_pulse: bool = True*) → None

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the pos\_pulse argument.

This function is blocking.

### Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **duration** – is the length of the impulse in seconds
- **pos\_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

**set\_t13\_socket** (*port: int, state: bool*) → None

Set the state of a SEV T13 power socket.

### Parameters

- **port** – is the socket number, one of *constants.T13\_SOCKET\_PORTS*
- **state** – is the desired on-state of the socket

**start()** → None

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

**stop()** → None

Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

**write** (*node\_id, value*) → None

Local wrapper for the OPC UA communication protocol write method.

### Parameters

- **node\_id** – the id of the node to read
- **value** – the value to write to the variable

```
class SupercubeConfiguration(namespace_index: int = 7)
Bases: object

Configuration dataclass for the Supercube devices.

clean_values()
    Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.

Parameters
    • fieldname – name of the field
    • value – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

namespace_index: int = 7
    Namespace of the OPC variables, typically this is 3 (coming from Siemens)

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

class SupercubeOpcUaCommunication(config)
Bases: hvl_ccb.comm.opc.OpcUaCommunication

Communication protocol specification for Supercube devices.

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

class SupercubeOpcUaCommunicationConfig(host: str, endpoint_name: str, port: int = 4845, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)
Bases: hvl_ccb.comm.opc.OpcUaCommunicationConfig

Communication protocol configuration for OPC UA, specifications for the Supercube devices.
```

**force\_value** (*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys** () → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults** () → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 4845**

Port of the OPC UA server to connect to.

**classmethod required\_keys** () → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**sub\_handler: hvl\_ccb.comm.opc.OpcUaSubHandler = <hvl\_ccb.dev.supercube2015.base.Super**

Subscription handler for data change events

**class SupercubeSubscriptionHandler**

Bases: *hvl\_ccb.comm.opc.OpcUaSubHandler*

OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.


**datachange\_notification** (*node: opcua.common.node.Node*, *val*, *data*)

In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.

**Parameters**

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

## hvl\_ccb.dev.supercube2015.constants

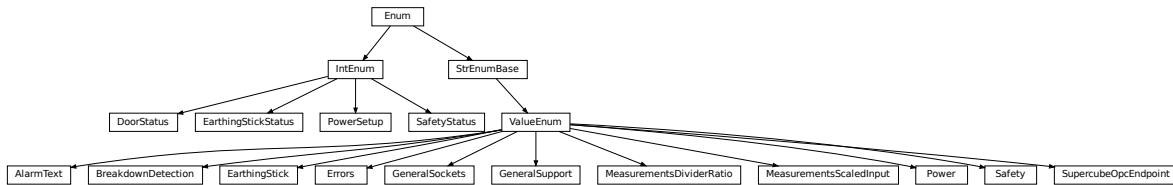

Constants, variable names for the Supercube OPC-connected devices.

```
class AlarmText (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the `AlarmText.get()` method to retrieve the enum of an alarm number.

```

Alarm0 = 'No Alarm.'

Alarm1 = 'STOP Safety switch 1 error'
Alarm10 = 'STOP Earthing stick 2 error'
Alarm11 = 'STOP Earthing stick 3 error'
Alarm12 = 'STOP Earthing stick 4 error'
Alarm13 = 'STOP Earthing stick 5 error'
Alarm14 = 'STOP Earthing stick 6 error'
Alarm17 = 'STOP Source switch error'
Alarm19 = 'STOP Fence 1 error'
Alarm2 = 'STOP Safety switch 2 error'
Alarm20 = 'STOP Fence 2 error'
Alarm21 = 'STOP Control error'
Alarm22 = 'STOP Power outage'
Alarm3 = 'STOP Emergency Stop 1'
Alarm4 = 'STOP Emergency Stop 2'
Alarm5 = 'STOP Emergency Stop 3'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error'
```

```
classmethod get(alarm: int)
```

Get the attribute of this enum for an alarm number.

**Parameters** `alarm` – the alarm number

**Returns** the enum for the desired alarm number

```
not_defined = 'NO ALARM TEXT DEFINED'

class BreakdownDetection(value=<object object>, names=None, module=None, type=None,
                           start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Node ID strings for the breakdown detection.

activated = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.Breakdowndetection.connect'
Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the system or not.

reset = 'hvl-ipc.WINAC.Support6OutA'
Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

triggered = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.Breakdowndetection.triggered'
Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using the safety circuit state, therefore no method is implemented to read this out directly.

class DoorStatus(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for doors.

closed = 2
Door is closed, but not locked.

error = 4
Door has an error or was opened in locked state (either with emergency stop or from the inside).

inactive = 0
not enabled in Supercube HMI setup, this door is not supervised.

locked = 3
Door is closed and locked (safe state).

open = 1
Door is open.

class EarthingStick(value=<object object>, names=None, module=None, type=None, start=1,
                     boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the earthing in manual mode.

classmethod manual(number: int)
Get the manual enum attribute for an earthing stick number.

Parameters number – the earthing stick (1..6)

Returns the manual enum

manual_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.MANUAL'
manual_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.MANUAL'
manual_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.MANUAL'
manual_4 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.MANUAL'
manual_5 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.MANUAL'
manual_6 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.MANUAL'
```

```

status_1_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1 CLOSE'
status_1_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1 CONNECT'
status_1_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1 OPEN'
status_2_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2 CLOSE'
status_2_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2 CONNECT'
status_2_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2 OPEN'
status_3_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3 CLOSE'
status_3_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3 CONNECT'
status_3_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3 OPEN'
status_4_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4 CLOSE'
status_4_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4 CONNECT'
status_4_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4 OPEN'
status_5_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5 CLOSE'
status_5_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5 CONNECT'
status_5_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5 OPEN'
status_6_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6 CLOSE'
status_6_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6 CONNECT'
status_6_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6 OPEN'

```

**classmethod status\_closed(number: int)**

Get the status enum attribute for an earthing stick number.

**Parameters** **number** – the earthing stick (1..6)

**Returns** the status enum

**classmethod status\_connected(number: int)**

Get the status enum attribute for an earthing stick number.

**Parameters** **number** – the earthing stick (1..6)

**Returns** the status enum

**classmethod status\_open(number: int)**

Get the status enum attribute for an earthing stick number.

**Parameters** **number** – the earthing stick (1..6)

**Returns** the status enum

```
class EarthingStickStatus(value=<object object>, names=None, module=None, type=None,
                         start=1, boundary=None)
```

Bases: aenum.IntEnum

Status of an earthing stick. These are the possible values in the status integer e.g. in EarthingStick. status\_1.

**closed = 1**

Earthing is closed (safe).

**error = 3**

Earthing is in error, e.g. when the stick did not close correctly or could not open.

```
inactive = 0
Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be
enabled in the Supercube HMI setup.

open = 2
Earthing is open (not safe).

class Errors(value=<object object>, names=None, module=None, type=None, start=1, bound-
ary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
Variable NodeID strings for information regarding error, warning and message handling.

quit = 'hvl-ipc.WINAC.SYSTEMSTATE.Faultconfirmation'
Writable boolean for the error quit button.

stop = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'
Boolean read-only variable telling if a stop is active.

stop_number = 'hvl-ipc.WINAC.SYSTEMSTATE.Errornumber'

class GeneralSockets(value=<object object>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
NodeID strings for the power sockets (3x T13 and 1xCEE16).

cee16 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.CEE16'
CEE16 socket (writeable boolean).

t13_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_1'
SEV T13 socket No. 1 (writable boolean).

t13_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_2'
SEV T13 socket No. 2 (writable boolean).

t13_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_3'
SEV T13 socket No. 3 (writable boolean).

class GeneralSupport(value=<object object>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
NodeID strings for the support inputs and outputs.

in_1_1 = 'hvl-ipc.WINAC.Support1InA'
in_1_2 = 'hvl-ipc.WINAC.Support1InB'
in_2_1 = 'hvl-ipc.WINAC.Support2InA'
in_2_2 = 'hvl-ipc.WINAC.Support2InB'
in_3_1 = 'hvl-ipc.WINAC.Support3InA'
in_3_2 = 'hvl-ipc.WINAC.Support3InB'
in_4_1 = 'hvl-ipc.WINAC.Support4InA'
in_4_2 = 'hvl-ipc.WINAC.Support4InB'
in_5_1 = 'hvl-ipc.WINAC.Support5InA'
in_5_2 = 'hvl-ipc.WINAC.Support5InB'
in_6_1 = 'hvl-ipc.WINAC.Support6InA'
```

```
in_6_2 = 'hvl-ipc.WINAC.Support6InB'
classmethod input (port, contact)
    Get the NodeID string for a support input.

    Parameters
        • port – the desired port (1..6)
        • contact – the desired contact at the port (1..2)

    Returns the node id string
```

```
out_1_1 = 'hvl-ipc.WINAC.Support1OutA'
out_1_2 = 'hvl-ipc.WINAC.Support1OutB'
out_2_1 = 'hvl-ipc.WINAC.Support2OutA'
out_2_2 = 'hvl-ipc.WINAC.Support2OutB'
out_3_1 = 'hvl-ipc.WINAC.Support3OutA'
out_3_2 = 'hvl-ipc.WINAC.Support3OutB'
out_4_1 = 'hvl-ipc.WINAC.Support4OutA'
out_4_2 = 'hvl-ipc.WINAC.Support4OutB'
out_5_1 = 'hvl-ipc.WINAC.Support5OutA'
out_5_2 = 'hvl-ipc.WINAC.Support5OutB'
out_6_1 = 'hvl-ipc.WINAC.Support6OutA'
out_6_2 = 'hvl-ipc.WINAC.Support6OutB'

classmethod output (port, contact)
    Get the NodeID string for a support output.
```

**Parameters**

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

**Returns** the node id string

```
class MeasurementsDividerRatio (value=<object object>, names=None, module=None,
                                type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

```
classmethod get (channel: int)
    Get the attribute for an input number.

    Parameters channel – the channel number (1..4)

    Returns the enum for the desired channel.
```

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.DivididerRatio'
class MeasurementsScaledInput (value=<object object>, names=None, module=None, type=None,
                                start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in [MeasurementsDividerRatio](#).

**classmethod** **get** (*channel: int*)

Get the attribute for an input number.

**Parameters** **channel** – the channel number (1..4)

**Returns** the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI1Volt'
input_2 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI2Volt'
input_3 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI3Volt'
input_4 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI4Volt'
```

**class** **Power** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

Variable NodeID strings concerning power data.

**current\_primary** = 'hvl-ipc.WINAC.SYSTEM\_INTERN.FUCurrentprim'

Primary current in ampere, measured by the frequency converter. (read-only)

**frequency** = 'hvl-ipc.WINAC.FU.Frequency'

Frequency converter output frequency. (read-only)

**setup** = 'hvl-ipc.WINAC.FU.TrafoSetup'

Power setup that is configured using the Supercube HMI. The value corresponds to the ones in [PowerSetup](#). (read-only)

**voltage\_max** = 'hvl-ipc.WINAC.FU.maxVoltagekV'

Maximum voltage allowed by the current experimental setup. (read-only)

**voltage\_primary** = 'hvl-ipc.WINAC.SYSTEM\_INTERN.FUVoltageprim'

Primary voltage in volts, measured by the frequency converter at its output. (read-only)

**voltage\_slope** = 'hvl-ipc.WINAC.FU.dUdt\_-1'

Voltage slope in V/s.

**voltage\_target** = 'hvl-ipc.WINAC.FU.SOLL'

Target voltage setpoint in V.

**class** **PowerSetup** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: [aenum.IntEnum](#)

Possible power setups corresponding to the value of variable [Power.setup](#).

**AC\_DoubleStage\_150kV** = 3

AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum voltage of 150kV.

**AC\_DoubleStage\_200kV** = 4

AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

**AC\_SingleStage\_100kV** = 2

AC voltage with MWB transformer set to 100kV maximum voltage.

**AC\_SingleStage\_50kV** = 1

AC voltage with MWB transformer set to 50kV maximum voltage.

**DC\_DoubleStage\_280kV = 7**

DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single stage transformer with Greinacher voltage doubling rectifier)

**DC\_SingleStage\_140kV = 6**

DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

**External = 0**

External power supply fed through blue CEE32 input using isolation transformer and safety switches of the Supercube, or using an external safety switch attached to the Supercube Type B.

**Internal = 5**

Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

```
class Safety(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

**horn = 'hvl-ipc.WINAC.SYSTEM\_INTERN.hornen'**

Writeable boolean to manually turn on or off the horn

**status\_error = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'****status\_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN'****status\_ready\_for\_red = 'hvl-ipc.WINAC.SYSTEMSTATE.ReadyForRed'**

Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in `SafetyStatus`.

**status\_red = 'hvl-ipc.WINAC.SYSTEMSTATE.RED'****switchto\_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN\_REQUEST'****switchto\_operate = 'hvl-ipc.WINAC.SYSTEMSTATE.switchon'**

Writable boolean for switching to Red Operate (locket, HV on) state.

**switchto\_ready = 'hvl-ipc.WINAC.SYSTEMSTATE.RED\_REQUEST'**

Writable boolean for switching to Red Ready (locked, HV off) state.

```
class SafetyStatus(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `aenum.IntEnum`

Safety status values that are possible states returned from `hvl_ccb.dev.supercube.base.Supercube.get_status()`. These values correspond to the states of the Supercube’s safety circuit statemachine.

**Error = 6**

System is in error mode.

**GreenNotReady = 1**

System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

**GreenReady = 2**

System is safe and all safety elements are in place to be able to switch to *ready*.

**Initializing = 0**

System is initializing or booting.

**QuickStop = 5**

Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

**RedOperate = 4**

System is locked in red state and in *operate* mode, i.e. high voltage on.

**RedReady = 3**

System is locked in red state and *ready* to go to *operate* mode.

```
class SupercubeOpcEndpoint (value=<object object>, names=None, module=None, type=None,
                             start=1, boundary=None)
```

Bases: [hvl\\_ccb.utils.enum.ValueEnum](#)

OPC Server Endpoint strings for the supercube variants.


A = 'OPC.SimaticNET.S7'

B = 'OPC.SimaticNET.S7'

**T13\_SOCKET\_PORTS = (1, 2, 3)**

Port numbers of SEV T13 power socket

## [hvl\\_ccb.dev.supercube2015.typ\\_a](#)

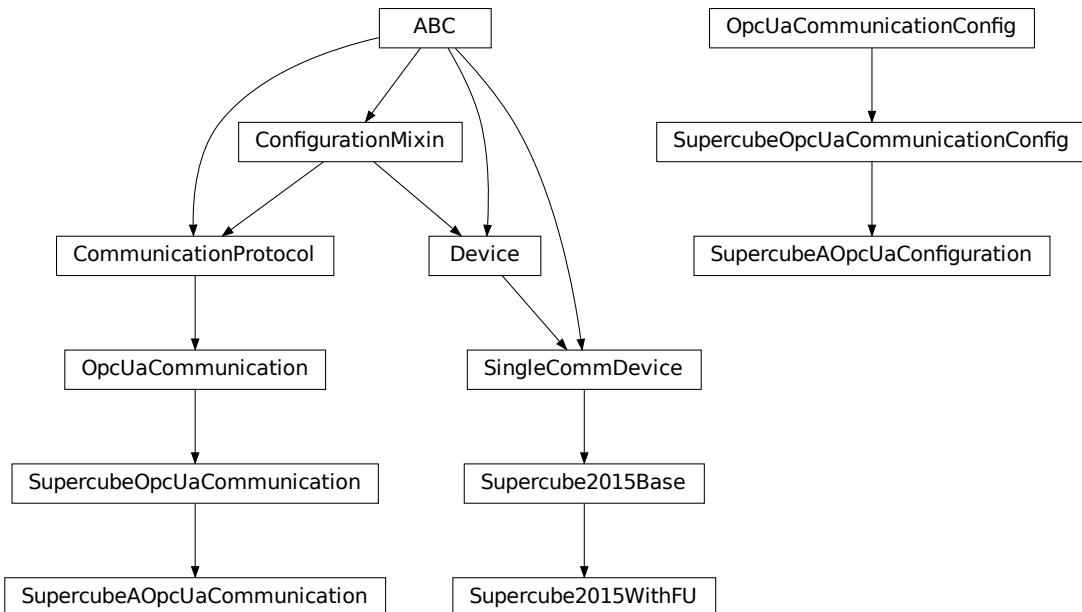

Supercube Typ A module.

```
class Supercube2015WithFU (com, dev_config=None)
```

Bases: [hvl\\_ccb.dev.supercube2015.base.Supercube2015Base](#)

Variant A of the Supercube with frequency converter.

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**fso\_reset()** → None

Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.

**get\_frequency()** → float

Read the electrical frequency of the current Supercube setup.

**Returns** the frequency in Hz

**get\_fso\_active()** → bool

Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.

**Returns** state of the FSO functionality

**get\_max\_voltage()** → float

Reads the maximum voltage of the setup and returns in V.

**Returns** the maximum voltage of the setup in V.

**get\_power\_setup()** → *hvl\_ccb.dev.supercube2015.constants.PowerSetup*

Return the power setup selected in the Supercube's settings.

**Returns** the power setup

**get\_primary\_current()** → float

Read the current primary current at the output of the frequency converter ( before transformer).

**Returns** primary current in A

**get\_primary\_voltage()** → float

Read the current primary voltage at the output of the frequency converter ( before transformer).

**Returns** primary voltage in V

**get\_target\_voltage()** → float

Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to *set\_target\_voltage()*.

**Returns** the setpoint voltage in V.

**set\_slope(*slope: float*)** → None

Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.

**Parameters** *slope* – voltage slope in V/s (0..15'000)

**set\_target\_voltage(*volt\_v: float*)** → None

Set the output voltage to a defined value in V.

**Parameters** *volt\_v* – the desired voltage in V

**class SupercubeAOpcUaCommunication(*config*)**

Bases: *hvl\_ccb.dev.supercube2015.base.SupercubeOpcUaCommunication*

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class SupercubeAOpcUaConfiguration(host: str, endpoint_name: str = "OPC.SimaticNET.S7", port: int = 4845, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object at 0x7f3720d9e4d0>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)
```

Bases: `hvl_ccb.dev.supercube2015.base.SuperCubeOpcUaCommunicationConfig`

**endpoint\_name:** `str = 'OPC.SimaticNET.S7'`  
Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

**force\_value** (`fieldname, value`)  
Forces a value to a dataclass field despite the class being frozen.  
NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


**Returns** a list of strings containing all required keys.

### Module contents

Supercube package with implementation for the old system version from 2015 based on Siemens WinAC soft-PLC on an industrial 32bit Windows computer.

## Submodules

### hvl\_ccb.dev.base

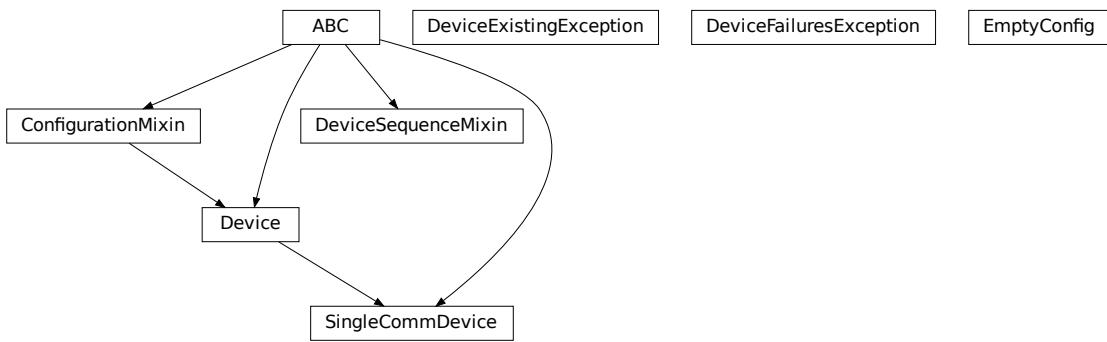

Module with base classes for devices.

**class Device(dev\_config=None)**

Bases: [hvl\\_ccb.configuration.ConfigurationMixin](#), [abc.ABC](#)

Base class for devices. Implement this class for a concrete device, such as measurement equipment or voltage sources.

Specifies the methods to implement for a device.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**abstract start()** → None

Start or restart this Device. To be implemented in the subclass.

**abstract stop()** → None

Stop this Device. To be implemented in the subclass.

**exception DeviceExistingException**

Bases: [Exception](#)

Exception to indicate that a device with that name already exists.

**exception DeviceFailuresException(failures: Dict[str, Exception], \*args)**

Bases: [Exception](#)

Exception to indicate that one or several devices failed.

**failures: Dict[str, Exception]**

A dictionary of named devices failures (exceptions).

**class DeviceSequenceMixin(devices: Dict[str, hvl\_ccb.dev.base.Device])**

Bases: [abc.ABC](#)

Mixin that can be used on a device or other classes to provide facilities for handling multiple devices in a sequence.

**add\_device** (*name: str, device: hvl\_ccb.dev.base.Device*) → None  
Add a new device to the device sequence.

**Parameters**

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

**Raises** *DeviceExistingException* –

**devices\_failed\_start:** Dict[str, hvl\_ccb.dev.base.Device]  
Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.  
Empty if *stop()* was called last; cf. *devices\_failed\_stop*.

**devices\_failed\_stop:** Dict[str, hvl\_ccb.dev.base.Device]  
Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.  
Empty if *start()* was called last; cf. *devices\_failed\_start*.

**get\_device** (*name: str*) → hvl\_ccb.dev.base.Device  
Get a device by name.

**Parameters** **name** – is the name of the device.

**Returns** the device object from this sequence.

**get\_devices** () → List[Tuple[str, hvl\_ccb.dev.base.Device]]  
Get list of name, device pairs according to current sequence.

**Returns** A list of tuples with name and device each.

**remove\_device** (*name: str*) → hvl\_ccb.dev.base.Device  
Remove a device from this sequence and return the device object.

**Parameters** **name** – is the name of the device.

**Returns** device object or *None* if such device was not in the sequence.

**Raises** **ValueError** – when device with given name was not found

**start** () → None  
Start all devices in this sequence in their added order.

**Raises** *DeviceFailuresException* – if one or several devices failed to start

**stop** () → None  
Stop all devices in this sequence in their reverse order.

**Raises** *DeviceFailuresException* – if one or several devices failed to stop

**class EmptyConfig**

Bases: object

Empty configuration dataclass that is the default configuration for a Device.

**clean\_values** ()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value** (*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

## Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys () → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults () → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys () → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**class SingleCommDevice (com, dev\_config=None)**

Bases: *hvl\_ccb.dev.base.Device*, abc.ABC

Base class for devices with a single communication protocol.

**property com**

Get the communication protocol of this device.

**Returns** an instance of CommunicationProtocol subtype

**abstract static default\_com\_cls () → Type[hvl\_ccb.comm.base.CommunicationProtocol]**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device


**start () → None**

Open the associated communication protocol.

**stop () → None**

Close the associated communication protocol.

## **hvl\_ccb.dev.crylas**

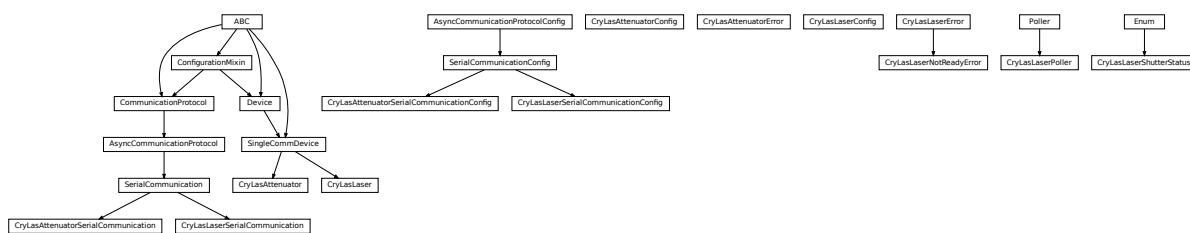

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: [https://www.crylas.de/products/pulsed\\_laser.html](https://www.crylas.de/products/pulsed_laser.html)

```
class CryLasAttenuator(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice
```

Device class for the CryLas laser attenuator.

```
property attenuation
```

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

```
set_attenuation(percent: Union[int, float]) → None
```

Set the percentage of attenuated light (inverse of set\_transmission). :param percent: percentage of attenuation, number between 0 and 100 :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

```
set_init_attenuation()
```

Sets the attenuation to its configured initial/default value

**Raises** `SerialCommunicationIOError` – when communication port is not opened

```
set_transmission(percent: Union[int, float]) → None
```

Set the percentage of transmitted light (inverse of set\_attenuation). :param percent: percentage of transmitted light :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

```
start() → None
```

Open the com, apply the config value ‘init\_attenuation’

**Raises** `SerialCommunicationIOError` – when communication port cannot be opened

```
property transmission
```

```
class CryLasAttenuatorConfig(init_attenuation: Union[int, float] = 0, response_sleep_time:
                                Union[int, float] = 1)
```

Bases: object

Device configuration dataclass for CryLas attenuator.

```
clean_values()
```

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

### Parameters

- **fieldname** – name of the field
- **value** – value to assign

```

init_attenuation: Union[int, float] = 0
is_configdataclass = True
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

response_sleep_time: Union[int, float] = 1

exception CryLasAttenuatorError
    Bases: Exception

    General error with the CryLas Attenuator.

class CryLasAttenuatorSerialCommunication(configuration)
    Bases: hvl\_ccb.comm.serial.SerialCommunication

    Specific communication protocol implementation for the CryLas attenuator. Already defines device-specific protocol parameters in config.

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

class CryLasAttenuatorSerialCommunicationConfig(terminator: bytes = b'', encoding: str = 'utf-8', encoding_error_handling: str = 'replace', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)
    Bases: hvl\_ccb.comm.serial.SerialCommunicationConfig

baudrate: int = 9600
    Baudrate for CryLas attenuator is 9600 baud

```

**bytesize:** Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8  
One byte is eight bits long

**force\_value** (fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'

CryLas attenuator does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1

CryLas attenuator uses one stop bit

**terminator:** bytes = b''

No terminator

**timeout:** Union[int, float] = 3

use 3 seconds timeout as default

**class CryLasLaser** (com, dev\_config=None)

Bases: hvl\_ccb.dev.base.SingleCommDevice

CryLas laser controller device class.

**class AnswersShutter** (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)

Bases: aenum.Enum

Standard answers of the CryLas laser controller to ‘Shutter’ command passed via *com*.

**CLOSED** = ‘Shutter inaktiv’

**OPENED** = ‘Shutter aktiv’

**class AnswersStatus** (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)

Bases: aenum.Enum

Standard answers of the CryLas laser controller to ‘STATUS’ command passed via *com*.

**ACTIVE** = ‘STATUS: Laser active’

```

HEAD = 'STATUS: Head ok'
INACTIVE = 'STATUS: Laser inactive'
READY = 'STATUS: System ready'
TEC1 = 'STATUS: TEC1 Regulation ok'
TEC2 = 'STATUS: TEC2 Regulation ok'

class LaserStatus(value=<object object>, names=None, module=None, type=None, start=1,
                  boundary=None)
Bases: aenum.Enum
Status of the CryLas laser
READY_ACTIVE = 2
READY_INACTIVE = 1
UNREADY_INACTIVE = 0
property is_inactive
property is_ready

class RepetitionRates(value=<object object>, names=None, module=None, type=None,
                      start=1, boundary=None)
Bases: aenum.IntEnum
Repetition rates for the internal software trigger in Hz
HARDWARE = 0
SOFTWARE_INTERNAL_SIXTY = 60
SOFTWARE_INTERNAL_TEN = 10
SOFTWARE_INTERNAL_TWENTY = 20

ShutterStatus
alias of hvl\_ccb.dev.crylas.CryLasLaserShutterStatus

close_shutter() → None
Close the laser shutter.

Raises
• SerialCommunicationIOError – when communication port is not opened
• CryLasLaserError – if success is not confirmed by the device

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_pulse_energy_and_rate() → Tuple[int, int]
Use the debug mode, return the measured pulse energy and rate.

Returns (energy in micro joule, rate in Hz)

Raises

```

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not answer the query

**laser\_off()** → None  
Turn the laser off.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**laser\_on()** → None  
Turn the laser on.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserNotReadyError* – if the laser is not ready to be turned on
- *CryLasLaserError* – if success is not confirmed by the device

**open\_shutter()** → None  
Open the laser shutter.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**set\_init\_shutter\_status()** → None  
Open or close the shutter, to match the configured shutter\_status.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**set\_pulse\_energy(energy: int)** → None  
Sets the energy of pulses (works only with external hardware trigger). Proceed with small energy steps, or the regulation may fail.

**Parameters** **energy** – energy in micro joule

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not confirm success

**set\_repetition\_rate(rate: Union[int, hvl\_ccb.dev.crylas.CryLasLaser.RepetitionRates])** → None  
Sets the repetition rate of the internal software trigger.

**Parameters** **rate** – frequency (Hz) as an integer

**Raises**

- *ValueError* – if rate is not an accepted value in RepetitionRates Enum
- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**start()** → None

Opens the communication protocol and configures the device.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened

**stop()** → None

Stops the device and closes the communication protocol.

**Raises**

- *SerialCommunicationIOError* – if com port is closed unexpectedly
- *CryLasLaserError* – if laser\_off() or close\_shutter() fail

**property target\_pulse\_energy**

**update\_laser\_status()** → None

Update the laser status to *LaserStatus.NOT\_READY* or *LaserStatus.INACTIVE* or *LaserStatus.ACTIVE*.

Note: laser never explicitly says that it is not ready (*LaserStatus.NOT\_READY*) in response to ‘STATUS’ command. It only says that it is ready (heated-up and implicitly inactive/off) or active (on). If it’s not either of these then the answer is *Answers.HEAD*. Moreover, the only time the laser explicitly says that its status is inactive (*Answers.INACTIVE*) is after issuing a ‘LASER OFF’ command.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**update\_repetition\_rate()** → None

Query the laser repetition rate.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**update\_shutter\_status()** → None

Update the shutter status (OPENED or CLOSED)

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**update\_target\_pulse\_energy()** → None

Query the laser pulse energy.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

**wait\_until\_ready()** → None

Block execution until the laser is ready

**Raises** *CryLasLaserError* – if the polling thread stops before the laser is ready

```
class CryLasLaserConfig(calibration_factor: Union[int, float] = 4.35, polling_period: Union[int, float] = 12, polling_timeout: Union[int, float] = 300, auto_laser_on: bool = True, init_shutter_status: Union[int, hvl_ccb.dev.crylas.CryLasLaserShutterStatus] = <CryLasLaserShutterStatus.CLOSED: 0>)
```

Bases: object

Device configuration dataclass for the CryLas laser controller.

**ShutterStatus**

alias of `hvl_ccb.dev.crylas.CryLasLaserShutterStatus`

**auto\_laser\_on:** bool = True

**calibration\_factor:** Union[int, float] = 4.35

**clean\_values()**

**force\_value** (fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**init\_shutter\_status:** Union[int, hvl\_ccb.dev.crylas.CryLasLaserShutterStatus] = 0

**is\_configdataclass** = True

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_period:** Union[int, float] = 12

**polling\_timeout:** Union[int, float] = 300

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception CryLasLaserError**

Bases: Exception

General error with the CryLas Laser.

**exception CryLasLaserNotReadyError**

Bases: `hvl_ccb.dev.crylas.CryLasLaserError`

Error when trying to turn on the CryLas Laser before it is ready.

**class CryLasLaserPoller** (spoll\_handler: Callable, check\_handler: Callable,  
check\_laser\_status\_handler: Callable, polling\_delay\_sec: Union[int,  
float] = 0, polling\_interval\_sec: Union[int, float] = 1, polling\_timeout\_sec:  
Optional[Union[int, float]] = None)

Bases: `hvl_ccb.dev.utils.Poller`

Poller class for polling the laser status until the laser is ready.

**Raises**

- `CryLasLaserError` – if the timeout is reached before the laser is ready

- *SerialCommunicationIOError* – when communication port is closed.

```
class CryLasLaserSerialCommunication(configuration)
    Bases: hvl_ccb.comm.serial.SerialCommunication
```

Specific communication protocol implementation for the CryLas laser controller. Already predefines device-specific protocol parameters in config.

```
READ_TEXT_SKIP_PREFIXES = ('>', 'MODE:')
```

Prefixes of lines that are skipped when read from the serial port.

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
query(cmd: str, prefix: str, post_cmd: Optional[str] = None) → str
```

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns the line in question.

#### Parameters

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer
- **post\_cmd** – optional additional command to send after the query

**Returns** line in question as a string

**Raises** *SerialCommunicationIOError* – when communication port is not opened

```
query_all(cmd: str, prefix: str)
```

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns a list of successive lines starting with prefix.

#### Parameters

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer

**Returns** line in question as a string

**Raises** *SerialCommunicationIOError* – when communication port is not opened

```
read() → str
```

Read first line of text from the serial port that does not start with any of *self.READ\_TEXT\_SKIP\_PREFIXES*.

**Returns** String read from the serial port; '' if there was nothing to read.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

```
class CryLasLaserSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str
= 'utf-8', encoding_error_handling: str =
'replace', wait_sec_read_text_nonempty:
Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int
= 10, port: Union[str, NoneType] = None, baudrate: int = 19200, parity: Union[str,
hvl_ccb.comm.serial.SerialCommunicationParity]
= <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
hvl_ccb.comm.serial.SerialCommunicationStopbits]
= <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
hvl_ccb.comm.serial.SerialCommunicationBytesize]
= <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int,
float] = 10)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate:** `int = 19200`

Baudrate for CryLas laser is 19200 baud

**bytesize:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8`

One byte is eight bits long

**force\_value** (`fieldname, value`)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** `Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'`

CryLas laser does not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1`

CryLas laser uses one stop bit

**terminator:** `bytes = b'\n'`

The terminator is LF

```

timeout: Union[int, float] = 10
    use 10 seconds timeout as default (a long timeout is needed!)


class CryLasLaserShutterStatus (value=<object object>, names=None, module=None,
    type=None, start=1, boundary=None)
Bases: aenum.Enum

Status of the CryLas laser shutter

CLOSED = 0
OPENED = 1

```

## [hvl\\_ccb.dev.ea\\_psi9000](#)

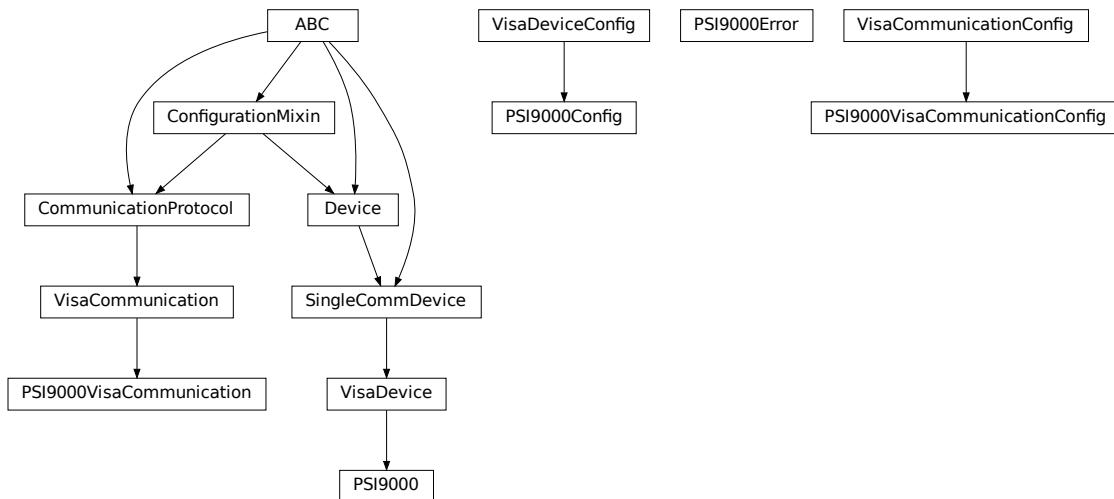

Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to know, all the testing was done with NI-Visa)

```

class PSI9000 (com: Union[hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunication,
    hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunicationConfig, dict], dev_config: Optional[Union[hvl_ccb.dev.ea_psi9000.PSI9000Config, dict]] = None)
Bases: hvl_ccb.dev.visa.VisaDevice

Elektro Automatik PSI 9000 power supply.

MS_NOMINAL_CURRENT = 2040
MS_NOMINAL_VOLTAGE = 80
SHUTDOWN_CURRENT_LIMIT = 0.1
SHUTDOWN_VOLTAGE_LIMIT = 0.1

check_master_slave_config() → None
    Checks if the master / slave configuration and initializes if successful

    Raises PSI9000Error – if master-slave configuration failed

```

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns** the VisaCommunication class

**get\_output() → bool**

Reads the current state of the DC output of the source. Returns True, if it is enabled, false otherwise.

**Returns** the state of the DC output

**get\_system\_lock() → bool**

Get the current lock state of the system. The lock state is true, if the remote control is active and false, if not.

**Returns** the current lock state of the device

**get\_ui\_lower\_limits() → Tuple[float, float]**

Get the lower voltage and current limits. A lower power limit does not exist.

**Returns** Umin in V, Imin in A

**get\_uip\_upper\_limits() → Tuple[float, float, float]**

Get the upper voltage, current and power limits.

**Returns** Umax in V, Imax in A, Pmax in W

**get\_voltage\_current\_setpoint() → Tuple[float, float]**

Get the voltage and current setpoint of the current source.

**Returns** Uset in V, Iset in A

**measure\_voltage\_current() → Tuple[float, float]**

Measure the DC output voltage and current

**Returns** Umeas in V, Imeas in A

**set\_lower\_limits(voltage\_limit: Optional[float] = None, current\_limit: Optional[float] = None)**

Set the lower limits for voltage and current. After writing the values a check is performed if the values are set correctly.

**Parameters**

- **voltage\_limit** – is the lower voltage limit in V
- **current\_limit** – is the lower current limit in A

**Raises** **PSI9000Error** – if the limits are out of range

**set\_output(target\_onstate: bool) → None**

Enables / disables the DC output.

**Parameters** **target\_onstate** – enable or disable the output power

**Raises** **PSI9000Error** – if operation was not successful

**set\_system\_lock(lock: bool) → None**

Lock / unlock the device, after locking the control is limited to this class unlocking only possible when voltage and current are below the defined limits

**Parameters** **lock** – True: locking, False: unlocking

---

**set\_upper\_limits** (*voltage\_limit: Optional[float] = None, current\_limit: Optional[float] = None, power\_limit: Optional[float] = None*) → None

Set the upper limits for voltage, current and power. After writing the values a check is performed if the values are set. If a parameter is left blank, the maximum configurable limit is set.

#### Parameters

- **voltage\_limit** – is the voltage limit in V
- **current\_limit** – is the current limit in A
- **power\_limit** – is the power limit in W

**Raises** `PSI9000Error` – if limits are out of range

**set\_voltage\_current** (*volt: float, current: float*) → None

Set voltage and current setpoints.

After setting voltage and current, a check is performed if writing was successful.

#### Parameters

- **volt** – is the setpoint voltage: 0..81.6 V (1.02 \* 0-80 V) (absolute max, can be smaller if limits are set)
- **current** – is the setpoint current: 0..2080.8 A (1.02 \* 0 - 2040 A) (absolute max, can be smaller if limits are set)

**Raises** `PSI9000Error` – if the desired setpoint is out of limits

**start** () → None

Start this device.

**stop** () → None

Stop this device. Turns off output and lock, if enabled.

**class PSI9000Config** (*spoll\_interval: Union[int, float] = 0.5, spoll\_start\_delay: Union[int, float] = 2, power\_limit: Union[int, float] = 43500, voltage\_lower\_limit: Union[int, float] = 0.0, voltage\_upper\_limit: Union[int, float] = 10.0, current\_lower\_limit: Union[int, float] = 0.0, current\_upper\_limit: Union[int, float] = 2040.0, wait\_sec\_system\_lock: Union[int, float] = 0.5, wait\_sec\_settings\_effect: Union[int, float] = 1, wait\_sec\_initialisation: Union[int, float] = 2*)

Bases: `hvl_ccb.dev.VisaDeviceConfig`

Elektro Automatik PSI 9000 power supply device class. The device is communicating over a VISA TCP socket.

Using this power supply, DC voltage and current can be supplied to a load with up to 2040 A and 80 V (using all four available units in parallel). The maximum power is limited by the grid, being at 43.5 kW available through the CEE63 power socket.

**clean\_values** () → None

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**current\_lower\_limit: Union[int, float] = 0.0**

Lower current limit in A, depending on the experimental setup.

**current\_upper\_limit: Union[int, float] = 2040.0**

Upper current limit in A, depending on the experimental setup.

**force\_value** (*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**power\_limit: Union[int, float] = 43500**

Power limit in W depending on the experimental setup. With 3x63A, this is 43.5kW. Do not change this value, if you do not know what you are doing. There is no lower power limit.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**voltage\_lower\_limit: Union[int, float] = 0.0**

Lower voltage limit in V, depending on the experimental setup.

**voltage\_upper\_limit: Union[int, float] = 10.0**

Upper voltage limit in V, depending on the experimental setup.

**wait\_sec\_initialisation: Union[int, float] = 2**

**wait\_sec\_settings\_effect: Union[int, float] = 1**

**wait\_sec\_system\_lock: Union[int, float] = 0.5**

**exception PSI9000Error**

Bases: Exception

Base error class regarding problems with the PSI 9000 supply.

**class PSI9000VisaCommunication(configuration)**

Bases: *hvl\_ccb.comm.visa.VisaCommunication*

Communication protocol used with the PSI 9000 power supply.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class PSI9000VisaCommunicationConfig(host: str, interface\_type: Union[str, *hvl\_ccb.comm.visa.VisaCommunicationConfig.InterfaceType*] = <InterfaceType.TCP/IP\_SOCKET: 1>, board: int = 0, port: int = 5025, timeout: int = 5000, chunk\_size: int = 204800, open\_timeout: int = 1000, write\_termination: str = '\n', read\_termination: str = '\n', visa\_backend: str = '')**

Bases: *hvl\_ccb.comm.visa.VisaCommunicationConfig*

Visa communication protocol config dataclass with specification for the PSI 9000 power supply.

**force\_value** (*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**interface\_type:** Union[str, *hvl\_ccb.comm.visa.VisaCommunicationConfig.InterfaceType*] = Interface type of the VISA connection, being one of `InterfaceType`.

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]


Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

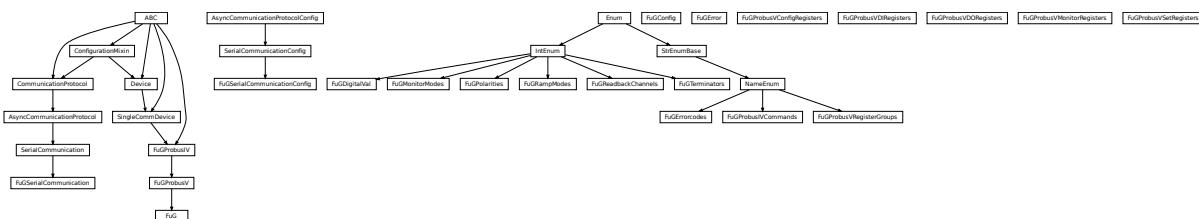**hvl\_ccb.dev.fug**

Device classes for “Probus V - ADDAT30” Interfaces which are used to control power supplies from FuG Elektronik GmbH

This interface is used for many FuG power units. Manufacturer homepage: <https://www.fug-elektronik.de>

The Professional Series of Power Supplies from FuG is a series of low, medium and high voltage direct current power supplies as well as capacitor chargers. The class `FuG` is tested with a HCK 800-20 000 in Standard Mode. The addressable mode is not implemented. Check the code carefully before using it with other devices. Manufacturer homepage: <https://www.fug-elektronik.de/netzgeraete/professional-series/>

The documentation of the interface from the manufacturer can be found here: [https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus\\_V.zip](https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus_V.zip)

The provided classes support the basic and some advanced commands. The commands for calibrating the power supplies are not implemented, as they are only for very special porpoises and should not be used by “normal” customers.

```
class FuG (com, dev_config=None)
Bases: hvl_ccb.dev.fug.FuGProbusV
```

FuG power supply device class.

The power supply is controlled over a FuG ADDA Interface with the PROBUS V protocol

**property config\_status**

Returns the registers for the registers with the configuration and status values

**Returns** FuGProbusVConfigRegisters

**property current**

Returns the registers for the current output

**Returns**

**property current\_monitor**

Returns the registers for the current monitor.

A typically usage will be “self.current\_monitor.value” to measure the output current

**Returns**

**property di**

Returns the registers for the digital inputs

**Returns** FuGProbusVDIRegisters

**identify\_device()** → None

Identify the device nominal voltage and current based on its model number.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**property max\_current**

Returns the maximal current which could provided within the test setup

**Returns**

**property max\_current\_hardware**

Returns the maximal current which could provided with the power supply

**Returns**

**property max\_voltage**

Returns the maximal voltage which could provided within the test setup

**Returns**

**property max\_voltage\_hardware**

Returns the maximal voltage which could provided with the power supply

**Returns**

**property on**

Returns the registers for the output switch to turn the output on or off

**Returns** FuGProbusVDORegisters

**property outx0**

Returns the registers for the digital output X0

**Returns** FuGProbusVDORegisters

**property outx1**

Returns the registers for the digital output X1

**Returns** FuGProbusVDORegisters

**property outx2**

Returns the registers for the digital output X2

**Returns** FuGProbusVDORegisters

**property outXCMD**

Returns the registers for the digital outputX-CMD

**Returns** FuGProbusVDORegisters

**start** (*max\_voltage=0, max\_current=0*) → None

Opens the communication protocol and configures the device.

**Parameters**

- **max\_voltage** – Configure here the maximal permissible voltage which is allowed in the given experimental setup
- **max\_current** – Configure here the maximal permissible current which is allowed in the given experimental setup

**property voltage**

Returns the registers for the voltage output

**Returns**

**property voltage\_monitor**

Returns the registers for the voltage monitor.

A typically usage will be “self.voltage\_monitor.value” to measure the output voltage

**Returns**

**class FuGConfig** (*wait\_sec\_stop\_commands: Union[int, float] = 0.5*)

Bases: object

Device configuration dataclass for FuG power supplies.

**clean\_values()**

**force\_value** (*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

wait_sec_stop_commands: Union[int, float] = 0.5
    Time to wait after subsequent commands during stop (in seconds)

class FuGDigitalVal(value)
    Bases: enum.IntEnum

    An enumeration.

    NO = 0
    OFF = 0
    ON = 1
    YES = 1

exception FuGError(*args, **kwargs)
    Bases: Exception

    Error with the FuG voltage source.

    errorcode: str
        Errorcode from the Probus, see documentation of Probus V chapter 5. Errors with three-digit errorcodes are thrown by this python module.

class FuGErrorcodes(value=<object object>, names=None, module=None, type=None, start=1,
                     boundary=None)
    Bases: hvl_ccb.utils.enum.NameEnum

The power supply can return an errorcode. These errorcodes are handled by this class. The original errorcodes from the source are with one or two digits, see documentation of Probus V chapter 5. All three-digit errorcodes are from this python module.

E0 = ('no error', 'standard response on each command')
E1 = ('no data available', 'Customer tried to read from GPIB but there were no data pr')
E10 = ('unknown SCPI command', 'This SCPI command is not implemented')
E100 = ('Command is not implemented', 'You tried to execute a command, which is not implem')
E106 = ('The rampstate is a read-only register', 'You tried to write data to the regis')
E11 = ('not allowed Trigger-on-Talk', 'Not allowed attempt to Trigger-on-Talk (~T1) wh')
E115 = ('The given index to select a digital value is out of range', 'Only integer val')
E12 = ('invalid argument in ~Tn command', 'Only ~T1 and ~T2 is implemented.')
E125 = ('The given index to select a ramp mode is out of range', 'Only integer values 1')
E13 = ('invalid N-value', 'Register > K8 contained an invalid value. Error code is out')
E135 = ('The given index to select the readback channel is out of range', 'Only integer')
E14 = ('register is write only', 'Some registers can only be writte to (i.e. > H0)')
E145 = ('The given value for the AD-conversion is unknown', 'Valid values for the ad-conv')
E15 = ('string too long', 'i.e. serial number string too long during calibration')
E155 = ('The given value to select a polarity is out range.', 'The value should be 0 or 1')


```

```

E16 = ('wrong checksum', 'checksum over command string was not correct, refer also to ...')
E165 = ('The given index to select the terminator string is out of range', '')
E2 = ('unknown register type', "No valid register type after '>'")
E206 = ('This status register is read-only', 'You tried to write data to this register')
E306 = ('The monitor register is read-only', 'You tried to write data to a monitor, whi... ')
E4 = ('invalid argument', 'The argument of the command was rejected .i.e. malformed num... ')
E5 = ('argument out of range', 'i.e. setvalue higher than type value')
E504 = ('Empty string as response', 'The connection is broken.')
E505 = ('The returned register is not the requested.', 'Maybe the connection is overbur... ')
E6 = ('register is read only', 'Some registers can only be read but not written to. (i... ')
E666 = ('You cannot overwrite the most recent error in the interface of the power supp... ')
E7 = ('Receive Overflow', 'Command string was longer than 50 characters.')
E8 = ('EEPROM is write protected', 'Write attempt to calibration data while the write p... ')
E9 = ('address error', 'A non addressed command was sent to ADDA while it was in address... ')
raise_()

class FuGMonitorModes(value)
Bases: enum.IntEnum

An enumeration.

T1MS = 1
    15 bit + sign, 1 ms integration time

T200MS = 6
    typ. 19 bit + sign, 200 ms integration time

T20MS = 3
    17 bit + sign, 20 ms integration time

T256US = 0
    14 bit + sign, 256 us integration time

T40MS = 4
    17 bit + sign, 40 ms integration time

T4MS = 2
    15 bit + sign, 4 ms integration time

T800MS = 7
    typ. 20 bit + sign, 800 ms integration time

T80MS = 5
    typ. 18 bit + sign, 80 ms integration time

class FuGPolarities(value)
Bases: enum.IntEnum

An enumeration.

NEGATIVE = 1
POSITIVE = 0

```

```
class FuGProbusIV(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice, abc.ABC
FuG Probus IV device class

Sends basic SCPI commands and reads the answer. Only the special commands and PROBUS IV instruction set is implemented.

command(command: hvl_ccb.dev.fug.FuGProbusIVCommands, value=None) → str

Parameters
    • command – one of the commands given within FuGProbusIVCommands
    • value – an optional value, depending on the command

Returns a String if a query was performed

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

output_off() → None
Switch DC voltage output off.

reset() → None
Reset of the interface: All setvalues are set to zero

abstract start()
Open the associated communication protocol.

stop() → None
Close the associated communication protocol.

class FuGProbusIVCommands(value=<object object>, names=None, module=None, type=None,
                           start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ADMODE = ('S', (<enum 'FuGMonitorModes'>, <class 'int'>))
CURRENT = ('I', (<class 'int'>, <class 'float'>))
EXECUTE = ('X', None)
EXECUTEONX = ('G', (<enum 'FuGDigitalVal'>, <class 'int'>))
    Wait for "X" to execute pending commands

ID = ('*IDN?', None)
OUTPUT = ('F', (<enum 'FuGDigitalVal'>, <class 'int'>))
POLARITY = ('P', (<enum 'FuGPolarities'>, <class 'int'>))
QUERY = ('?', None)
READBACKCHANNEL = ('N', (<enum 'FuGReadbackChannels'>, <class 'int'>))
RESET = ('=', None)
TERMINATOR = ('Y', (<enum 'FuGTerminators'>, <class 'int'>))
```

```

VOLTAGE = ('U', (<class 'int'>, <class 'float'>))
XOUTPUTS = ('R', <class 'int'>
    TODO: the possible values are limited to 0..13

class FuGProbusV(com, dev_config=None)
Bases: hvl_ccb.dev.fug.FuGProbusIV

FuG Probus V class which uses register based commands to control the power supplies

get_register(register: str) → str
    get the value from a register

        Parameters register – the register from which the value is requested

        Returns the value of the register as a String

set_register(register: str, value: Union[int, float, str]) → None
    generic method to set value to register

        Parameters
            • register – the name of the register to set the value
            • value – which should be written to the register

class FuGProbusVConfigRegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Configuration and Status values, acc. 4.2.5

property execute_on_x
    status of Execute-on-X

        Returns FuGDigitalVal of the status

property most_recent_error
    Reads the Error-Code of the most recent command

        Return FuGError
        Raises FuGError – if code is not “E0”

property readback_data
    Preselection of readout data for Trigger-on-Talk

        Returns index for the readback channel

property srq_mask
    SRQ-Mask, Service-Request Enable status bits for SRQ 0: no SRQ Bit 2: SRQ on change of status to CC
    Bit 1: SRQ on change to CV

        Returns representative integer value

property srq_status
    SRQ-Statusbyte output as a decimal number: Bit 2: PS is in CC mode Bit 1: PS is in CV mode

        Returns representative string

property status
    Statusbyte as a string of 0/1. Combined status (compatibel to Probus IV), MSB first: Bit 7: I-REG Bit
    6: V-REG Bit 5: ON-Status Bit 4: 3-Reg Bit 3: X-Stat (polarity) Bit 2: Cal-Mode Bit 1: unused Bit 0:
    SEL-D

        Returns string of 0/1

```

**property terminator**

Terminator character for answer strings from ADDA

**Returns** FuGTerminators

**class FuGProbusVDIRegisters** (*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Digital Inputs acc. 4.2.4

**property analog\_control**

**Returns** shows 1 if power supply is controlled by the analog interface

**property calibration\_mode**

**Returns** shows 1 if power supply is in calibration mode

**property cc\_mode**

**Returns** shows 1 if power supply is in CC mode

**property cv\_mode**

**Returns** shows 1 if power supply is in CV mode

**property digital\_control**

**Returns** shows 1 if power supply is digitally controlled

**property on**

**Returns** shows 1 if power supply ON

**property reg\_3**

For special applications.

**Returns** input from bit 3-REG

**property x\_stat**

**Returns** polarity of HVPS with polarity reversal

**class FuGProbusVDORegisters** (*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Digital outputs acc. 4.2.2

**property out**

Status of the output according to the last setting. This can differ from the actual state if output should only pulse.

**Returns** FuGDigitalVal

**property status**

Returns the actual value of output. This can differ from the set value if pulse function is used.

**Returns** FuGDigitalVal

**class FuGProbusVMonitorRegisters** (*fug, super\_register: hvl\_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Analog monitors acc. 4.2.3

**property adc\_mode**

The programmed resolution and integration time of the AD converter

**Returns** FuGMonitorModes

---

**property value**  
Value from the monitor.

**Returns** a float value in V or A

**property value\_raw**  
uncalibrated raw value from AD converter

**Returns** float value from ADC

```
class FuGProbusVRegisterGroups (value=<object object>, names=None, module=None,
                                 type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
```

An enumeration.

```
CONFIG = 'K'
INPUT = 'D'
MONITOR_I = 'M1'
MONITOR_V = 'M0'
OUTPUTONCMD = 'BON'
OUTPUTX0 = 'B0'
OUTPUTX1 = 'B1'
OUTPUTX2 = 'B2'
OUTPUTXCMD = 'BX'
SETCURRENT = 'S1'
SETVOLTAGE = 'S0'
```

```
class FuGProbusVSetRegisters (fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object
```

Setvalue control acc. 4.2.1 for the voltage and the current output

**property actualsetvalue**  
The actual valid set value, which depends on the ramp function.

**Returns** actual valid set value

**property high\_resolution**  
Status of the high resolution mode of the output.

**Return 0** normal operation  
**Return 1** High Res. Mode

**property rampmode**  
The set ramp mode to control the setvalue.

**Returns** the mode of the ramp as instance of FuGRampModes

**property ramprate**  
The set ramp rate in V/s.

**Returns** ramp rate in V/s

**property rampstate**  
Status of ramp function.

**Return 0** if final setvalue is reached

**Return 1** if still ramping up

**property setvalue**

For the voltage or current output this setvalue was programmed.

**Returns** the programmed setvalue

**class FuGRampModes (value)**

Bases: enum.IntEnum

An enumeration.

**FOLLOWRAMP = 1**

Follow the ramp up- and downwards

**IMMEDIATELY = 0**

Standard mode: no ramp

**ONLYUPWARDSOFFTOZERO = 4**

Follow the ramp up- and downwards, if output is OFF set value is zero

**RAMPUPWARDS = 2**

Follow the ramp only upwards, downwards immediately

**SPECIALRAMPUPWARDS = 3**

Follow a special ramp function only upwards

**class FuGReadbackChannels (value)**

Bases: enum.IntEnum

An enumeration.

**CURRENT = 1**

**FIRMWARE = 5**

**RATEDCURRENT = 4**

**RATEDVOLTAGE = 3**

**SN = 6**

**STATUSBYTE = 2**

**VOLTAGE = 0**

**class FuGSerialCommunication (configuration)**

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for FuG power supplies. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query (command: str) → str**

Send a command to the interface and handle the status message. Eventually raises an exception.

**Parameters** **command** – Command to send

**Raises** **FuGError** – if the connection is broken or the error from the power source itself

**Returns** Answer from the interface or empty string

```
class FuGSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-
8', encoding_error_handling: str = 'replace',
wait_sec_read_text_nonempty: Union[int, float] =
0.5, default_n_attempts_read_text_nonempty:
int = 10, port: Union[str, NoneType] = None,
baudrate: int = 9600, parity: Union[str,
hvl_ccb.comm.serial.SerialCommunicationParity]
= <SerialCommunicationPar-
ity.NONE: 'N'>, stopbits: Union[int,
hvl_ccb.comm.serial.SerialCommunicationStopbits]
= <SerialCommunicationStop-
bits.ONE: 1>, bytesize: Union[int,
hvl_ccb.comm.serial.SerialCommunicationBytesize] =
<SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
Union[int, float] = 3)
Bases: hvl_ccb.comm.serial.SerialCommunicationConfig
```

**baudrate: int = 9600**  
Baudrate for FuG power supplies is 9600 baud

**bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8**  
One byte is eight bits long

**default\_n\_attempts\_read\_text\_nonempty: int = 10**  
default number of attempts to read a non-empty text

**force\_value (fieldname, value)**  
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys () → Sequence[str]**  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults () → Dict[str, object]**  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'**  
FuG does not use parity

**classmethod required\_keys () → Sequence[str]**  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1**  
FuG uses one stop bit

**terminator: bytes = b'\n'**  
The terminator is LF

```

timeout: Union[int, float] = 3
    use 3 seconds timeout as default

wait_sec_read_text_nonempty: Union[int, float] = 0.5
    default time to wait between attempts of reading a non-empty text


class FuGTerminators(value)
    Bases: enum.IntEnum

    An enumeration.

    CR = 3
    CRLF = 0
    LF = 2
    LFCR = 1

```

## **hvl\_ccb.dev.heinzinger**

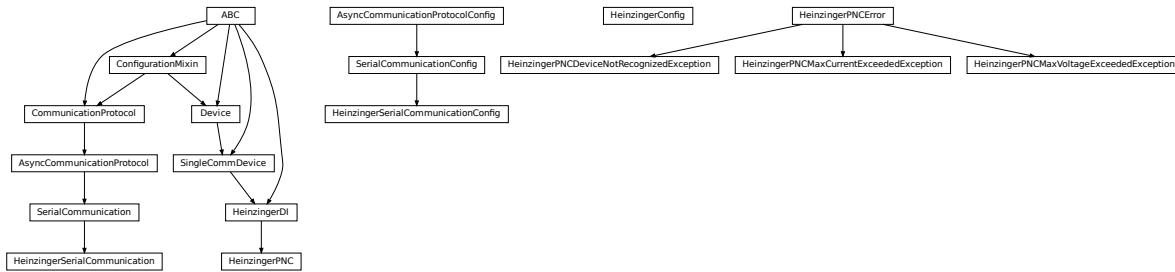

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Manufacturer homepage: <https://www.heinzinger.com/products/accessories-and-more/digital-interfaces/>

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class **HeinzingerPNC** is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCCap. Manufacturer homepage: <https://www.heinzinger.com/products/high-voltage/universal-high-voltage-power-supplies/>

```

class HeinzingerConfig(default_number_of_recordings:
                           Union[int,
                                 hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum]] = 1, num-
                           ber_of_decimals: int = 6, wait_sec_stop_commands: Union[int, float] =
                           0.5)

```

Bases: object

Device configuration dataclass for Heinzinger power supplies.

```
class RecordingsEnum(value)
```

Bases: enum.IntEnum

An enumeration.

```
EIGHT = 8
```

```
FOUR = 4
```

```
ONE = 1
```

```
SIXTEEN = 16
```

```

TWO = 2

clean_values()
default_number_of_recordings: Union[int, hvl_ccb.dev.heinzinger.HeinzingerConfig.Recording]
force_value(fieldname, value)
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define post_force_value method with same signature as this method to do extra processing
after value has been forced on fieldname.

Parameters

- fieldname – name of the field
- value – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

number_of_decimals: int = 6

classmethod optional_defaults() → Dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
instantiation.

Returns a list of strings containing all required keys.

wait_sec_stop_commands: Union[int, float] = 0.5
Time to wait after subsequent commands during stop (in seconds)

class HeinzingerDI(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice, abc.ABC
Heinzinger Digital Interface I/II device class
Sends basic SCPI commands and reads the answer. Only the standard instruction set from the manual is implemented.

class OutputStatus(value)
Bases: enum.IntEnum
Status of the voltage output
OFF = 0
ON = 1
UNKNOWN = -1

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

```

```
static default_com_cls()
Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

get_current() → float
Queries the set current of the Heinzinger PNC (not the measured current!).

    Raises SerialCommunicationIOError – when communication port is not opened

get_interface_version() → str
Queries the version number of the digital interface.

    Raises SerialCommunicationIOError – when communication port is not opened

get_number_of_recordings() → int
Queries the number of recordings the device is using for average value calculation.

    Returns int number of recordings

    Raises SerialCommunicationIOError – when communication port is not opened

get_serial_number() → str
Ask the device for its serial number and returns the answer as a string.

    Returns string containing the device serial number

    Raises SerialCommunicationIOError – when communication port is not opened

get_voltage() → float
Queries the set voltage of the Heinzinger PNC (not the measured voltage!).

    Raises SerialCommunicationIOError – when communication port is not opened

measure_current() → float
Ask the Device to measure its output current and return the measurement result.

    Returns measured current as float

    Raises SerialCommunicationIOError – when communication port is not opened

measure_voltage() → float
Ask the Device to measure its output voltage and return the measurement result.

    Returns measured voltage as float

    Raises SerialCommunicationIOError – when communication port is not opened

output_off() → None
Switch DC voltage output off and updates the output status.

    Raises SerialCommunicationIOError – when communication port is not opened

output_on() → None
Switch DC voltage output on and updates the output status.

    Raises SerialCommunicationIOError – when communication port is not opened

property output_status
reset_interface() → None
Reset of the digital interface; only Digital Interface I: Power supply is switched to the Local-Mode (Manual operation)

    Raises SerialCommunicationIOError – when communication port is not opened
```

---

**set\_current** (*value: Union[int, float]*) → None  
Sets the output current of the Heinzinger PNC to the given value.

**Parameters** **value** – current expressed in *self.unit\_current*

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_number\_of\_recordings** (*value: Union[int, hvl\_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum]*) → None  
Sets the number of recordings the device is using for average value calculation. The possible values are 1, 2, 4, 8 and 16.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_voltage** (*value: Union[int, float]*) → None  
Sets the output voltage of the Heinzinger PNC to the given value.

**Parameters** **value** – voltage expressed in *self.unit\_voltage*

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**abstract start()**  
Opens the communication protocol.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened.

**stop()** → None  
Stop the device. Closes also the communication protocol.

**class HeinzingerPNC** (*com, dev\_config=None*)  
Bases: *hvl\_ccb.dev.heinzinger.HeinzingerDI*

Heinzinger PNC power supply device class.

The power supply is controlled over a Heinzinger Digital Interface I/II

**class UnitCurrent** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)  
Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

An enumeration.

**A = 3**

**UNKNOWN = 1**

**mA = 2**

**class UnitVoltage** (*value=<object object>, names=None, module=None, type=None, start=1, boundary=None*)  
Bases: *hvl\_ccb.utils.enum.AutoNumberNameEnum*

An enumeration.

**UNKNOWN = 1**

**v = 2**

**kV = 3**

**identify\_device()** → None  
Identify the device nominal voltage and current based on its serial number.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**property max\_current**

**property max\_current\_hardware**

```
property max_voltage
property max_voltage_hardware
set_current (value: Union[int, float]) → None
    Sets the output current of the Heinzinger PNC to the given value.

    Parameters value – current expressed in self.unit_current
    Raises SerialCommunicationIOError – when communication port is not opened

set_voltage (value: Union[int, float]) → None
    Sets the output voltage of the Heinzinger PNC to the given value.

    Parameters value – voltage expressed in self.unit_voltage
    Raises SerialCommunicationIOError – when communication port is not opened

start () → None
    Opens the communication protocol and configures the device.

property unit_current
property unit_voltage

exception HeinzingerPNCDeviceNotRecognizedException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError
    Error indicating that the serial number of the device is not recognized.

exception HeinzingerPNCError
    Bases: Exception
    General error with the Heinzinger PNC voltage source.

exception HeinzingerPNCMaxCurrentExceededException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError
    Error indicating that program attempted to set the current to a value exceeding ‘max_current’.

exception HeinzingerPNCMaxVoltageExceededException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError
    Error indicating that program attempted to set the voltage to a value exceeding ‘max_voltage’.

class HeinzingerSerialCommunication (configuration)
    Bases: hvl_ccb.comm.serial.SerialCommunication
    Specific communication protocol implementation for Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

static config_cls ()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class
```

```
class HeinzingerSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str
= 'utf-8', encoding_error_handling: str =
'replace', wait_sec_read_text_nonempty:
Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int
= 40, port: Union[str, NoneType] = None,
baudrate: int = 9600, parity: Union[str,
hvl_ccb.comm.serial.SerialCommunicationParity]
= <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
hvl_ccb.comm.serial.SerialCommunicationStopbits]
= <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
hvl_ccb.comm.serial.SerialCommunicationBytesize]
= <SerialCommunicationBytesize.EIGHTBITS:
8>, timeout: Union[int, float] = 3)
Bases: hvl_ccb.comm.serial.SerialCommunicationConfig
```

**baudrate: int = 9600**

Baudrate for Heinzinger power supplies is 9600 baud

**bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8**

One byte is eight bits long

**default\_n\_attempts\_read\_text\_nonempty: int = 40**

increased to 40 default number of attempts to read a non-empty text

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field

- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'**

Heinzinger does not use parity

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1**


Heinzinger uses one stop bit

```
terminator: bytes = b'\n'
The terminator is LF

timeout: Union[int, float] = 3
use 3 seconds timeout as default

wait_sec_read_text_nonempty: Union[int, float] = 0.5
default time to wait between attempts of reading a non-empty text
```

## hvl\_ccb.dev.labjack

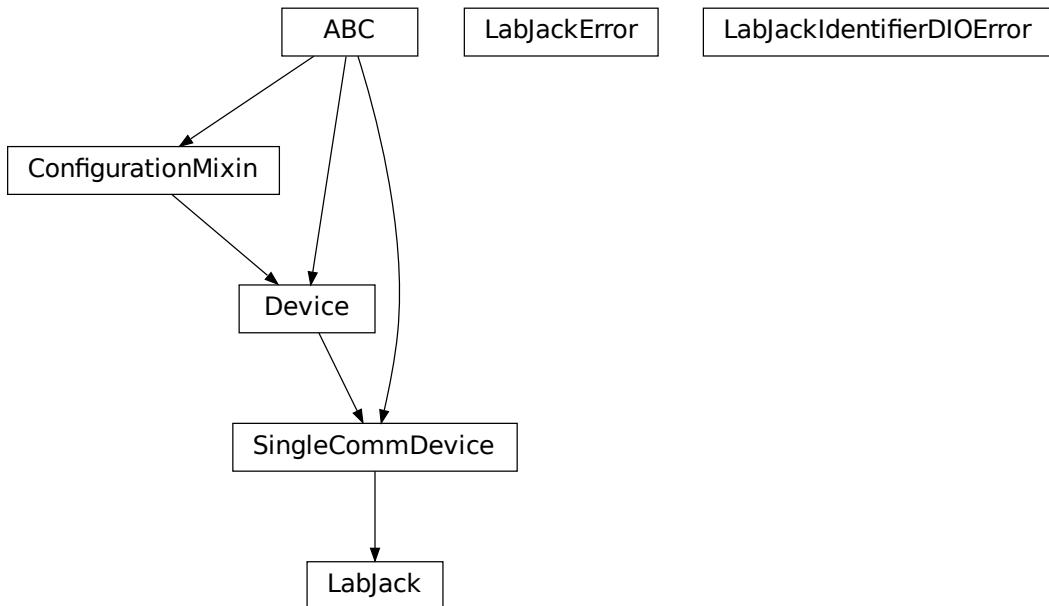

Labjack Device for hvl\_ccb. Originally developed and tested for LabJack T7-PRO.

Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

```
class LabJack(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice
```

LabJack Device.

This class is tested with a LabJack T7-Pro and should also work with T4 and T7 devices communicating through the LJM Library. Other or older hardware versions and variants of LabJack devices are not supported.

```
class AInRange(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.StrEnumBase
```

An enumeration.

```
ONE = 1.0
```

```

ONE_HUNDREDTH = 0.01
ONE_TENTH = 0.1
TEN = 10.0
property value

class CalMicroAmpere(value=<object object>, names=None, module=None, type=None,
                           start=1, boundary=None)
    Bases: aenum.Enum
    Pre-defined microampere (uA) values for calibration current source query.

TEN = '10uA'
TWO_HUNDRED = '200uA'

class CjcType(value=<object object>, names=None, module=None, type=None, start=1, bound-
ary=None)
    Bases: hvl_ccb.utils.enum.NameEnum
    CJC slope and offset
    internal = (1, 0)
    lm34 = (55.56, 255.37)

DIOChannel
    alias of hvl_ccb._dev.labjack.TSeriesDIOChannel

class DIOStatus(value=<object object>, names=None, module=None, type=None, start=1,
                      boundary=None)
    Bases: aenum.IntEnum
    State of a digital I/O channel.

HIGH = 1
LOW = 0

class DeviceType(value=<object object>, names=None, module=None, type=None, start=1,
                           boundary=None)
    Bases: hvl_ccb.utils.enum.AutoNumberNameEnum
    LabJack device types.

    Can be also looked up by ambiguous Product ID (p_id) or by instance name: ``python
    LabJackDeviceType(4) is LabJackDeviceType('T4') ``

ANY = 1
T4 = 2
T7 = 3
T7_PRO = 4

classmethod get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
                                         List[hvl_ccb._dev.labjack.DeviceType]]
    Get LabJack device type instance via LabJack product ID.

    Note: Product ID is not unambiguous for LabJack devices.

    Parameters p_id – Product ID of a LabJack device
    Returns Instance or list of instances of LabJackDeviceType
    Raises ValueError – when Product ID is unknown

```

```

class TemperatureUnit(value=<object object>, names=None, module=None, type=None,
                      start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.NameEnum

Temperature unit (to be returned)

C = 1
F = 2
K = 0

class ThermocoupleType(value=<object object>, names=None, module=None, type=None,
                           start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.NameEnum

Thermocouple type; NONE means disable thermocouple mode.

C = 30
E = 20
J = 21
K = 22
NONE = 0
PT100 = 40
PT1000 = 42
PT500 = 41
R = 23
S = 25
T = 24

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_ain(*channels: int) → Union[float, Sequence[float]]
Read currently measured value (voltage, resistance, ...) from one or more of analog inputs.

Parameters channels – AIN number or numbers (0..254)

Returns the read value (voltage, resistance, ...) as float or 'tuple of them in case multiple channels given

get_cal_current_source(name: Union[str, CalMicroAmpere]) → float
This function will return the calibration of the chosen current source, this ist not a measurement!

The value was stored during fabrication.

Parameters name – ‘200uA’ or ‘10uA’ current source

Returns calibration of the chosen current source in ampere

get_digital_input(address: Union[str, hvl\_ccb.\_dev.labjack.TSeriesDIOChannel]) →
                    hvl\_ccb.dev.labjack.LabJack.DIOStatus
Get the value of a digital input.

```

allowed names for T7 (Pro): FIO0 - FIO7, EIO0 - EIO 7, CIO0- CIO3, MIO0 - MIO2 :param address: name of the output -> ‘FIO0’ :return: HIGH when *address* DIO is high, and LOW when *address* DIO is low

**get\_product\_id()** → int

This function returns the product ID reported by the connected device.

Attention: returns 7 for both T7 and T7-Pro devices!

**Returns** integer product ID of the device

**get\_product\_name(*force\_query\_id=False*)** → str

This function will return the product name based on product ID reported by the device.

Attention: returns “T7” for both T7 and T7-Pro devices!

**Parameters** **force\_query\_id** – boolean flag to force *get\_product\_id* query to device instead of using cached device type from previous queries.

**Returns** device name string, compatible with *LabJack.DeviceType*

**get\_product\_type(*force\_query\_id: bool = False*)** → hvl\_ccb.\_dev.labjack.DeviceType

This function will return the device type based on reported device type and in case of unambiguity based on configuration of device’s communication protocol (e.g. for “T7” and “T7\_PRO” devices), or, if not available first matching.

**Parameters** **force\_query\_id** – boolean flag to force *get\_product\_id* query to device instead of using cached device type from previous queries.

**Returns** *DeviceType* instance

**Raises** *LabJackIdentifierDIOError* – when read Product ID is unknown

**get\_sbus\_rh(*number: int*)** → float

Read the relative humidity value from a serial SBUS sensor.

**Parameters** **number** – port number (0..22)

**Returns** relative humidity in %RH

**get\_sbus\_temp(*number: int*)** → float

Read the temperature value from a serial SBUS sensor.

**Parameters** **number** – port number (0..22)

**Returns** temperature in Kelvin

**get\_serial\_number()** → int

Returns the serial number of the connected LabJack.

**Returns** Serial number.

**read\_resistance(*channel: int*)** → float

Read resistance from specified channel.

**Parameters** **channel** – channel with resistor

**Returns** resistance value with 2 decimal places

**read\_thermocouple(*pos\_channel: int*)** → float

Read the temperature of a connected thermocouple.

**Parameters** **pos\_channel** – is the AIN number of the positive pin

**Returns** temperature in specified unit

**set\_ain\_differential** (*pos\_channel*: int, *differential*: bool) → None

Sets an analog input to differential mode or not. T7-specific: For base differential channels, positive must be even channel from 0-12 and negative must be positive+1. For extended channels 16-127, see Mux80 datasheet.

#### Parameters

- **pos\_channel** – is the AIN number (0..12)
- **differential** – True or False

**Raises** *LabJackError* – if parameters are unsupported

**set\_ain\_range** (*channel*: int, *vrange*: Union[Real, AInRange]) → None

Set the range of an analog input port.

#### Parameters

- **channel** – is the AIN number (0..254)
- **vrange** – is the voltage range to be set

**set\_ain\_resistance** (*channel*: int, *vrange*: Union[Real, AInRange], *resolution*: int) → None

Set the specified channel to resistance mode. It utilized the 200uA current source of the LabJack.

#### Parameters

- **channel** – channel that should measure the resistance
- **vrange** – voltage range of the channel
- **resolution** – resolution index of the channel T4: 0-5, T7: 0-8, T7-Pro 0-12

**set\_ain\_resolution** (*channel*: int, *resolution*: int) → None

Set the resolution index of an analog input port.

#### Parameters

- **channel** – is the AIN number (0..254)
- **resolution** – is the resolution index within 0...`get\_product\_type().ain\_max\_resolution` range; 0 will set the resolution index to default value.

**set\_ain\_thermocouple** (*pos\_channel*: int, *thermocouple*: Union[None, str, ThermocoupleType],  
                          *cjc\_address*: int = 60050, *cjc\_type*: Union[str, CjcType] = <*Cjc-*  
                          *Type.internal*: (1, 0)>, *vrange*: Union[Real, AInRange] = <*AIn-*  
                          *Range.ONE\_HUNDREDTH*: '0.01'>, *resolution*: int = 10, *unit*: Union[str,  
                          *TemperatureUnit*] = <*TemperatureUnit.K*: 0>) → None

Set the analog input channel to thermocouple mode.

#### Parameters

- **pos\_channel** – is the analog input channel of the positive part of the differential pair
- **thermocouple** – None to disable thermocouple mode, or string specifying the thermocouple type
- **cjc\_address** – modbus register address to read the CJC temperature
- **cjc\_type** – determines cjc slope and offset, ‘internal’ or ‘lm34’
- **vrange** – measurement voltage range
- **resolution** – resolution index (T7-Pro: 0-12)
- **unit** – is the temperature unit to be returned (‘K’, ‘C’ or ‘F’)

**Raises** *LabJackError* – if parameters are unsupported

**set\_digital\_output** (*address: str, state: Union[int, DIOStatus]*) → None  
Set the value of a digital output.

#### Parameters

- **address** – name of the output -> ‘FIO0’
- **state** – state of the output -> *DIOStatus* instance or corresponding *int* value

**start()** → None  
Start the Device.

**stop()** → None  
Stop the Device.

#### exception LabJackError

Bases: Exception


Errors of the LabJack device.

#### exception LabJackIdentifierDIOError

Bases: Exception

Error indicating a wrong DIO identifier

### hvl\_ccb.dev.mbw973

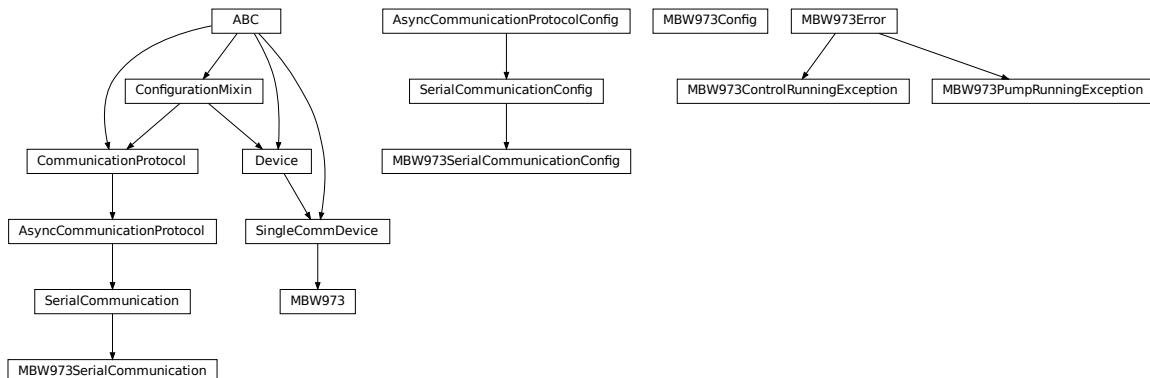

Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

**class MBW973 (com, dev\_config=None)**  
Bases: *hvl\_ccb.dev.base.SingleCommDevice*

MBW 973 dew point mirror device class.

**static config\_cls()**  
Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**  
Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**is\_done()** → bool

Poll status of the dew point mirror and return True, if all measurements are done.

**Returns** True, if all measurements are done; False otherwise.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**read**(*cast\_type*: Type = <class 'str'>)

Read value from *self.com* and cast to *cast\_type*. Raises *ValueError* if read text (*str*) is not convertible to *cast\_type*, e.g. to *float* or to *int*.

**Returns** Read value of *cast\_type* type.

**read\_float()** → float

Convenience wrapper for *self.read()*, with typing hint for return value.

**Returns** Read *float* value.

**read\_int()** → int

Convenience wrapper for *self.read()*, with typing hint for return value.

**Returns** Read *int* value.

**read\_measurements()** → Dict[str, float]

Read out measurement values and return them as a dictionary.

**Returns** Dictionary with values.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**set\_measuring\_options**(*humidity*: bool = True, *sf6\_purity*: bool = False) → None

Send measuring options to the dew point mirror.

#### Parameters

- **humidity** – Perform humidity test or not?
- **sf6\_purity** – Perform SF6 purity test or not?

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**start()** → None

Start this device. Opens the communication protocol and retrieves the set measurement options from the device.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened.

**start\_control()** → None

Start dew point control to acquire a new value set.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**stop()** → None

Stop the device. Closes also the communication protocol.

**write**(*value*) → None

Send *value* to *self.com*.

**Parameters** **value** – Value to send, converted to *str*.

**Raises** *SerialCommunicationIOError* – when communication port is not opened

**class MBW973Config**(*polling\_interval*: Union[int, float] = 2)

Bases: object

Device configuration dataclass for MBW973.

**clean\_values ()**

**force\_value (fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys () → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults () → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_interval: Union[int, float] = 2**

Polling period for *is\_done* status queries [in seconds].

**classmethod required\_keys () → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception MBW973ControlRunningException**

Bases: *hvl\_ccb.dev.mbw973.MBW973Error*

Error indicating there is still a measurement running, and a new one cannot be started.

**exception MBW973Error**

Bases: *Exception*

General error with the MBW973 dew point mirror device.

**exception MBW973PumpRunningException**

Bases: *hvl\_ccb.dev.mbw973.MBW973Error*

Error indicating the pump of the dew point mirror is still recovering gas, unable to start a new measurement.

**class MBW973SerialCommunication (configuration)**

Bases: *hvl\_ccb.comm.serial.SerialCommunication*

Specific communication protocol implementation for the MBW973 dew point mirror. Already predefines device-specific protocol parameters in config.

**static config\_cls ()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class MBW973SerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8', encoding_error_handling: str = 'replace', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)
Bases: hvl_ccb.comm.serial.SerialCommunicationConfig
```

**baudrate: int = 9600**  
Baudrate for MBW973 is 9600 baud

**bytesize: Union[int, hvl\_ccb.comm.serial.SerialCommunicationBytesize] = 8**  
One byte is eight bits long

**force\_value(fieldname, value)**  
Forces a value to a dataclass field despite the class being frozen.  
NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity: Union[str, hvl\_ccb.comm.serial.SerialCommunicationParity] = 'N'**  
MBW973 does not use parity

**classmethod required\_keys() → Sequence[str]**  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


**Returns** a list of strings containing all required keys.

**stopbits: Union[int, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1**  
MBW973 does use one stop bit

**terminator: bytes = b'\r'**  
The terminator is only CR

**timeout: Union[int, float] = 3**  
use 3 seconds timeout as default

## hvl\_ccb.dev.newport

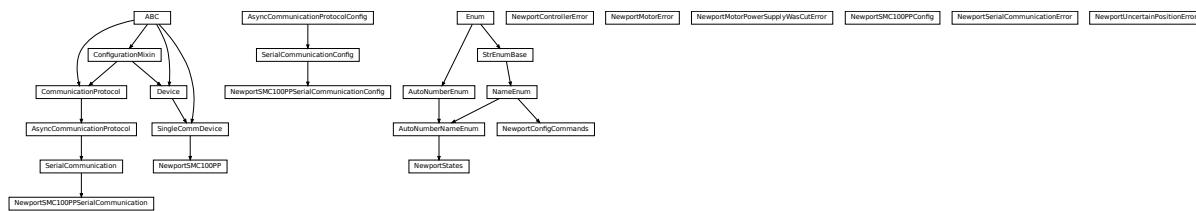

Device class for Newport SMC100PP stepper motor controller with serial communication.

The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/f/smci00-single-axis-dc-or-stepper-motion-controller>

```
class NewportConfigCommands (value=<object object>, names=None, module=None, type=None,
                           start=1, boundary=None)
```

Bases: `hvl_ccb.utils.enum.NameEnum`

Commands predefined by the communication protocol of the SMC100PP

```
AC = 'acceleration'
BA = 'backlash_compensation'
BH = 'hysteresis_compensation'
FRM = 'micro_step_per_full_step_factor'
FRS = 'motion_distance_per_full_step'
HT = 'home_search_type'
JR = 'jerk_time'
OH = 'home_search_velocity'
OT = 'home_search_timeout'
QIL = 'peak_output_current_limit'
SA = 'rs485_address'
SL = 'negative_software_limit'
SR = 'positive_software_limit'
VA = 'velocity'
VB = 'base_velocity'
ZX = 'stage_configuration'
```

```
exception NewportControllerError
```

Bases: `Exception`

Error with the Newport controller.

```
exception NewportMotorError
```

Bases: `Exception`

Error with the Newport motor.

```
exception NewportMotorPowerSupplyWasCutError
Bases: Exception
```

Error with the Newport motor after the power supply was cut and then restored, without interrupting the communication with the controller.

```
class NewportSMC100PP (com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice
```

Device class of the Newport motor controller SMC100PP

```
class MotorErrors (value=<object object>, names=None, module=None, type=None, start=1,
boundary=None)
```

Bases: aenum.Enum

Possible motor errors reported by the motor during get\_state().

```
DC_VOLTAGE_TOO_LOW = 3
FOLLOWING_ERROR = 6
HOMING_TIMEOUT = 5
NED_END_OF_TURN = 11
OUTPUT_POWER_EXCEEDED = 2
PEAK_CURRENT_LIMIT = 9
POS_END_OF_TURN = 10
RMS_CURRENT_LIMIT = 8
SHORT_CIRCUIT = 7
WRONG_ESP_STAGE = 4
```

```
class StateMessages (value=<object object>, names=None, module=None, type=None, start=1,
boundary=None)
```

Bases: aenum.Enum

Possible messages returned by the controller on get\_state() query.

```
CONFIG = '14'
DISABLE_FROM_JOGGING = '3E'
DISABLE_FROM_MOVING = '3D'
DISABLE_FROM_READY = '3C'
HOMING_FROM_RS232 = '1E'
HOMING_FROM_SMC = '1F'
JOGGING_FROM_DISABLE = '47'
JOGGING_FROM_READY = '46'
MOVING = '28'
NO_REF_ESP_STAGE_ERROR = '10'
NO_REF_FROM_CONFIG = '0C'
NO_REF_FROM_DISABLED = '0D'
NO_REF_FROM_HOMING = '0B'
NO_REF_FROM_JOGGING = '11'
```

```

NO_REF_FROM_MOVING = '0F'
NO_REF_FROM_READY = '0E'
NO_REF_FROM_RESET = '0A'
READY_FROM_DISABLE = '34'
READY_FROM_HOMING = '32'
READY_FROM_JOGGING = '35'
READY_FROM_MOVING = '33'

```

**States**

alias of `hvl_ccb.dev.newport.NewportStates`

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**exit\_configuration (add: Optional[int] = None) → None**

Exit the CONFIGURATION state and go back to the NOT REFERENCED state. All configuration parameters are saved to the device's memory.

**Parameters** `add` – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**get\_acceleration (add: Optional[int] = None) → Union[int, float]**

Leave the configuration state. The configuration parameters are saved to the device's memory.

**Parameters** `add` – controller address (1 to 31)

**Returns** acceleration (preset units/s^2), value between 1e-6 and 1e12

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**get\_controller\_information (add: Optional[int] = None) → str**

Get information on the controller name and driver version

**Parameters** `add` – controller address (1 to 31)

**Returns** controller information

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained

- *NewportControllerError* – if the controller reports an error

**get\_motor\_configuration** (*add: Optional[int] = None*) → Dict[str, float]

Query the motor configuration and returns it in a dictionary.

**Parameters** **add** – controller address (1 to 31)

**Returns** dictionary containing the motor's configuration

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_move\_duration** (*dist: Union[int, float], add: Optional[int] = None*) → float

Estimate the time necessary to move the motor of the specified distance.

**Parameters**

- **dist** – distance to travel
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_negative\_software\_limit** (*add: Optional[int] = None*) → Union[int, float]

Get the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

**Parameters** **add** – controller address (1 to 31)

**Returns** negative software limit (preset units), value between -1e12 and 0

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**get\_position** (*add: Optional[int] = None*) → float

Returns the value of the current position.

**Parameters** **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportUncertainPositionError* – if the position is ambiguous

**get\_positive\_software\_limit** (*add: Optional[int] = None*) → Union[int, float]

Get the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

**Parameters** `add` – controller address (1 to 31)

**Returns** positive software limit (preset units), value between 0 and 1e12

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**get\_state** (`add: int = None`) → `StateMessages`

Check on the motor errors and the controller state

**Parameters** `add` – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error
- `NewportMotorError` – if the motor reports an error

**Returns** state message from the device (member of `StateMessages`)

**go\_home** (`add: Optional[int] = None`) → None

Move the motor to its home position.

**Parameters** `add` – controller address (1 to 31), defaults to self.address

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**go\_to\_configuration** (`add: Optional[int] = None`) → None

This method is executed during start(). It can also be executed after a reset(). The controller is put in CONFIG state, where configuration parameters can be changed.

**Parameters** `add` – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**initialize** (`add: Optional[int] = None`) → None

Puts the controller from the NOT\_REF state to the READY state. Sends the motor to its “home” position.

**Parameters** `add` – controller address (1 to 31)

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**move\_to\_absolute\_position** (*pos: Union[int, float]*, *add: Optional[int] = None*) → None  
Move the motor to the specified position.

**Parameters**

- **pos** – target absolute position (affected by the configured offset)
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**move\_to\_relative\_position** (*pos: Union[int, float]*, *add: Optional[int] = None*) → None  
Move the motor of the specified distance.

**Parameters**

- **pos** – distance to travel (the sign gives the direction)
- **add** – controller address (1 to 31), defaults to self.address

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**reset** (*add: Optional[int] = None*) → None

Resets the controller, equivalent to a power-up. This puts the controller back to NOT REFERENCED state, which is necessary for configuring the controller.

**Parameters** **add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**set\_acceleration** (*acc: Union[int, float]*, *add: Optional[int] = None*) → None

Leave the configuration state. The configuration parameters are saved to the device's memory.

**Parameters**

- **acc** – acceleration (preset units/s<sup>2</sup>), value between 1e-6 and 1e12
- **add** – controller address (1 to 31)

**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

**set\_motor\_configuration** (*add: Optional[int] = None*, *config: Optional[dict] = None*) → None  
Set the motor configuration. The motor must be in CONFIG state.

**Parameters**

- **add** – controller address (1 to 31)
- **config** – dictionary containing the motor's configuration

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**set\_negative\_software\_limit** (*lim: Union[int, float]*, *add: Optional[int] = None*) → None

Set the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

**Parameters**

- **lim** – negative software limit (preset units), value between -1e12 and 0
- **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**set\_positive\_software\_limit** (*lim: Union[int, float]*, *add: Optional[int] = None*) → None

Set the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

**Parameters**

- **lim** – positive software limit (preset units), value between 0 and 1e12
- **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**start()**

Opens the communication protocol and applies the config.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened

**stop()** → None

Stop the device. Close the communication protocol.

**stop\_motion** (*add: Optional[int] = None*) → None

Stop a move in progress by decelerating the positioner immediately with the configured acceleration until it stops. If a controller address is provided, stops a move in progress on this controller, else stops the moves on all controllers.

**Parameters** **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained

- *NewportControllerError* – if the controller reports an error

**wait\_until\_motor\_initialized**(*add*: *Optional[int]* = *None*) → *None*

Wait until the motor leaves the HOMING state (at which point it should have arrived to the home position).

**Parameters** *add* – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

```
class NewportSMC100PPConfig(address: int = 1, user_position_offset: Union[int,
    float] = 23.987, screw_scaling: Union[int, float] = 1,
    exit_configuration_wait_sec: Union[int, float] = 5, move_wait_sec:
    Union[int, float] = 1, acceleration: Union[int, float] = 10, backlash_compensation:
    Union[int, float] = 0, hysteresis_compensation:
    Union[int, float] = 0.015, micro_step_per_full_step_factor:
    int = 100, motion_distance_per_full_step: Union[int,
    float] = 0.01, home_search_type: Union[int,
    hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] =
    <HomeSearch.HomeSwitch: 2>, jerk_time: Union[int, float] = 0.04,
    home_search_velocity: Union[int, float] = 4, home_search_timeout:
    Union[int, float] = 27.5, home_search_polling_interval:
    Union[int, float] = 1, peak_output_current_limit: Union[int,
    float] = 0.4, rs485_address: int = 2, negative_software_limit:
    Union[int, float] = -23.5, positive_software_limit: Union[int,
    float] = 25, velocity: Union[int, float] = 4, base_velocity:
    Union[int, float] = 0, stage_configuration: Union[int,
    hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageConfig] =
    <EspStageConfig.EnableEspStageCheck: 3>)
```

Bases: *object*

Configuration dataclass for the Newport motor controller SMC100PP.

```
class EspStageConfig(value=<object object>, names=None, module=None, type=None,
    start=1, boundary=None)
```

Bases: *aenum.IntEnum*

Different configurations to check or not the motor configuration upon power-up.

**DisableEspStageCheck** = 1

**EnableEspStageCheck** = 3

**UpdateEspStageInfo** = 2

```
class HomeSearch(value=<object object>, names=None, module=None, type=None, start=1,
    boundary=None)
```

Bases: *aenum.IntEnum*

Different methods for the motor to search its home position during initialization.

**CurrentPosition** = 1

**EndOfRunSwitch** = 4

**EndOfRunSwitch\_and\_Index** = 3

**HomeSwitch** = 2

```

HomeSwitch_and_Index = 0
acceleration: Union[int, float] = 10
address: int = 1
backlash_compensation: Union[int, float] = 0
base_velocity: Union[int, float] = 0
clean_values()
exit_configuration_wait_sec: Union[int, float] = 5
force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

```

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
home_search_polling_interval: Union[int, float] = 1
```

```
home_search_timeout: Union[int, float] = 27.5
```

```
home_search_type: Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] = 1
```

```
home_search_velocity: Union[int, float] = 4
```

```
hysteresis_compensation: Union[int, float] = 0.015
```

```
is_configdataclass = True
```

```
jerk_time: Union[int, float] = 0.04
```

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

```
micro_step_per_full_step_factor: int = 100
```

```
motion_distance_per_full_step: Union[int, float] = 0.01
```

```
property motor_config
```

Gather the configuration parameters of the motor into a dictionary.

**Returns** dict containing the configuration parameters of the motor

```
move_wait_sec: Union[int, float] = 1
```

```
negative_software_limit: Union[int, float] = -23.5
```

```
classmethod optional_defaults() → Dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```
peak_output_current_limit: Union[int, float] = 0.4
```

```
positive_software_limit: Union[int, float] = 25
```

```
post_force_value(fieldname, value)
```

```
classmethod required_keys() → Sequence[str]
```

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

```
rs485_address: int = 2
```

```
screw_scaling: Union[int, float] = 1
```

```
stage_configuration: Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageCon
```

```
user_position_offset: Union[int, float] = 23.987
```

```
velocity: Union[int, float] = 4
```

```
class NewportSMC100PPSerialCommunication(configuration)
```

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

```
class ControllerErrors(value=<object object>, names=None, module=None, type=None,  
                      start=1, boundary=None)
```

Bases: [aenum.Enum](#)

Possible controller errors with values as returned by the device in response to sent commands.

```
ADDR_INCORRECT = 'B'
```

```
CMD_EXEC_ERROR = 'V'
```

```
CMD_NOT_ALLOWED = 'D'
```

```
CMD_NOT_ALLOWED_CC = 'X'
```

```
CMD_NOT_ALLOWED_CONFIGURATION = 'I'
```

```
CMD_NOT_ALLOWED_DISABLE = 'J'
```

```
CMD_NOT_ALLOWED_HOMING = 'L'
```

```
CMD_NOT_ALLOWED_MOVING = 'M'
```

```
CMD_NOT_ALLOWED_NOT_REFERENCED = 'H'
```

```
CMD_NOT_ALLOWED_PP = 'W'
```

```
CMD_NOT_ALLOWED_READY = 'K'
```

```
CODE_OR_ADDR_INVALID = 'A'
```

```
COM_TIMEOUT = 'S'
```

```
DISPLACEMENT_OUT_OF_LIMIT = 'G'
```

```
EEPROM_ACCESS_ERROR = 'U'
```

```
ESP_STAGE_NAME_INVALID = 'F'
```

```
HOME_STARTED = 'E'
```

```
NO_ERROR = '@'
```

```
PARAM_MISSING_OR_INVALID = 'C'
```

```
POSITION_OUT_OF_LIMIT = 'N'
```

**check\_for\_error** (*add: int*) → None

Ask the Newport controller for the last error it recorded.

This method is called after every command or query.

**Parameters** **add** – controller address (1 to 31)

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query** (*add: int, cmd: str, param: Optional[Union[int, float, str]] = None*) → str

Send a query to the controller, read the answer, and check for errors. The prefix add+cmd is removed from the answer.

**Parameters**

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

**Returns** the answer from the device without the prefix

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**query\_multiple** (*add: int, cmd: str, prefixes: List[str]*) → List[str]

Send a query to the controller, read the answers, and check for errors. The prefixes are removed from the answers.

**Parameters**

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **prefixes** – prefixes of each line expected in the answer

**Returns** list of answers from the device without prefix

**Raises**

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

**read\_text()** → str

Read one line of text from the serial port, and check for presence of a null char which indicates that the motor power supply was cut and then restored. The input buffer may hold additional data afterwards, since only one line is read.

This method uses `self.access_lock` to ensure thread-safety.

**Returns** String read from the serial port; ‘’ if there was nothing to read.

**Raises**

- `SerialCommunicationIOError` – when communication port is not opened
- `NewportMotorPowerSupplyWasCutError` – if a null char is read

**send\_command** (`add: int, cmd: str, param: Optional[Union[int, float, str]] = None`) → None

Send a command to the controller, and check for errors.

**Parameters**

- `add` – the controller address (1 to 31)
- `cmd` – the command to be sent
- `param` – optional parameter (int/float/str) appended to the command

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

**send\_stop** (`add: int`) → None

Send the general stop ST command to the controller, and check for errors.

**Parameters** `add` – the controller address (1 to 31)

**Returns** ControllerErrors reported by Newport Controller

**Raises**

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained

```
class NewportSMC100PPSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding_error_handling: str = 'replace', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 57600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 10)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate: int = 57600**

Baudrate for Heinzinger power supplies is 9600 baud

**bytesize:** Union[int, [hvl\\_ccb.comm.serial.SerialCommunicationBytesize](#)] = 8  
One byte is eight bits long

**force\_value** (fieldname, value)  
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys** () → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults** () → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** Union[str, [hvl\\_ccb.comm.serial.SerialCommunicationParity](#)] = 'N'  
Heinzinger does not use parity

**classmethod required\_keys** () → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**stopbits:** Union[int, [hvl\\_ccb.comm.serial.SerialCommunicationStopbits](#)] = 1  
Heinzinger uses one stop bit

**terminator:** bytes = b'\r\n'  
The terminator is CR/LF

**timeout:** Union[int, float] = 10  
use 10 seconds timeout as default

#### exception NewportSerialCommunicationError

Bases: Exception

Communication error with the Newport controller.

**class NewportStates** (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: [hvl\\_ccb.utils.enum.AutoNumberNameEnum](#)

States of the Newport controller. Certain commands are allowed only in certain states.

**CONFIG** = 3

**DISABLE** = 6

**HOMING** = 2


**JOGGING** = 7

**MOVING** = 5

**NO\_REF** = 1

```
READY = 4
exception NewportUncertainPositionError
    Bases: Exception
    Error with the position of the Newport motor.
```

## hvl\_ccb.dev.pfeiffer\_tpg

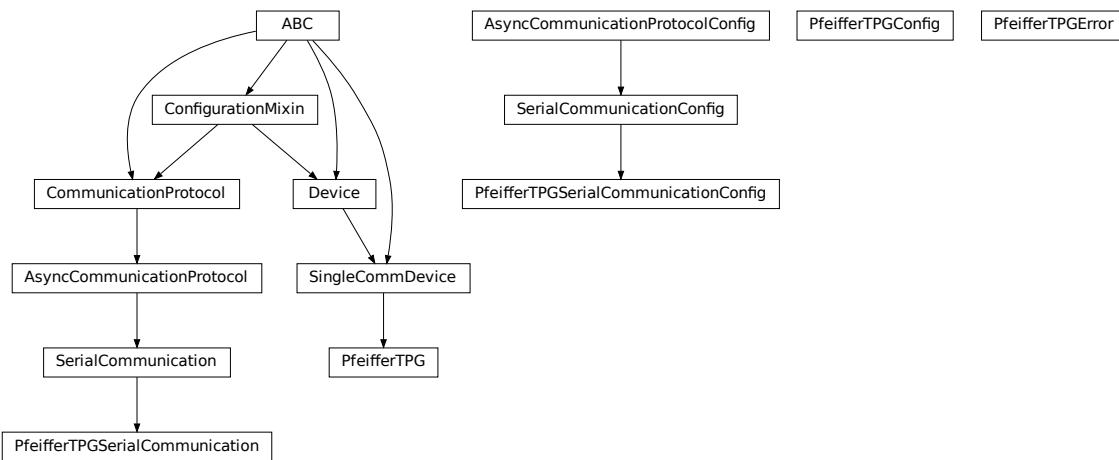

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/> measurement/activeline/controllers/

```
class PfeifferTPG (com, dev_config=None)
    Bases: hvl_ccb.dev.base.SingleCommDevice
```

Pfeiffer TPG control unit device class

```
class PressureUnits (value=<object object>, names=None, module=None, type=None, start=1,
                     boundary=None)
    Bases: hvl_ccb.utils.enum.NameEnum
```

Enum of available pressure units for the digital display. “0” corresponds either to bar or to mbar depending on the TPG model. In case of doubt, the unit is visible on the digital display.

```
Micron = 3
Pascal = 2
Torr = 1
Volt = 5
bar = 0
hPascal = 4
mbar = 0
```

---

```
class SensorStatus(value)
```

Bases: enum.IntEnum

An enumeration.

```
Identification_error = 6
```

```
No_sensor = 5
```

```
Ok = 0
```

```
Overrange = 2
```

```
Sensor_error = 3
```

```
Sensor_off = 4
```

```
Underrange = 1
```

```
class SensorTypes(value)
```

Bases: enum.Enum

An enumeration.

```
CMR = 4
```

```
IKR = 2
```

```
IKR11 = 2
```

```
IKR9 = 2
```

```
IMR = 5
```

```
None = 7
```

```
PBR = 6
```

```
PKR = 3
```

```
TPR = 1
```

```
noSENSOR = 7
```

```
noSen = 7
```

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
static default_com_cls()
```

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

```
get_full_scale_mbar() → List[Union[int, float]]
```

Get the full scale range of the attached sensors

**Returns** full scale range values in mbar, like [0.01, 1, 0.1, 1000, 50000, 10]

**Raises**

- **SerialCommunicationIOError** – when communication port is not opened

- **PfeifferTPGError** – if command fails

`get_full_scale_unitless() → List[int]`

Get the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

**Returns** list of full scale range values, like [0, 1, 3, 3, 2, 0]

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

`identify_sensors() → None`

Send identification request TID to sensors on all channels.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

`measure(channel: int) → Tuple[str, float]`

Get the status and measurement of one sensor

**Parameters** `channel` – int channel on which the sensor is connected, with  $1 \leq \text{channel} \leq \text{number\_of\_sensors}$

**Returns** measured value as float if measurement successful, sensor status as string if not

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

`measure_all() → List[Tuple[str, float]]`

Get the status and measurement of all sensors (this command is not available on all models)

**Returns** list of measured values as float if measurements successful, and or sensor status as strings if not

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

`property number_of_sensors`

`set_display_unit(unit: Union[str, hvl_ccb.dev.pfeiffer_tpg.PfeifferTPG.PressureUnits]) → None`

Set the unit in which the measurements are shown on the display.

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

`set_full_scale_mbar(fsr: List[Union[int, float]]) → None`

Set the full scale range of the attached sensors (in unit mbar)

**Parameters** `fsr` – full scale range values in mbar, for example [0.01, 1000]

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**set\_full\_scale\_unitless** (*fsr: List[int]*) → None  
Set the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

**Parameters** *fsr* – list of full scale range values, like [0, 1, 3, 3, 2, 0]

**Raises**

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

**start()** → None  
Start this device. Opens the communication protocol, and identify the sensors.

**Raises** *SerialCommunicationIOError* – when communication port cannot be opened

**stop()** → None  
Stop the device. Closes also the communication protocol.

**property unit**  
The pressure unit of readings is always mbar, regardless of the display unit.

```
class PfeifferTPGConfig(model: Union[str, hvl_ccb.dev.pfeiffer_tpg.PfeifferTPGConfig.Model] = <Model.TPG25xA: {1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7, 0.1: 8}>)
```

Bases: *object*  
Device configuration dataclass for Pfeiffer TPG controllers.

```
class Model(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: *hvl\_ccb.utils.enum.NameEnum*  
An enumeration.

```
TPG25xA = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6}
```

```
TPGx6x = {0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8}
```

```
is_valid_scale_range_reversed_str(v: str) → bool
```

Check if given string represents a valid reversed scale range of a model.

**Parameters** *v* – Reversed scale range string.  
**Returns** *True* if valid, *False* otherwise.

**clean\_values()**

**force\_value** (*fieldname, value*)  
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]  
Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

```
model: Union[str, hvl_ccb.dev.pfeiffer_tpg.PfeifferTPGConfig.Model] = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7, 0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8}
```

**classmethod optional\_defaults()** → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception PfeifferTPGError**

Bases: Exception

Error with the Pfeiffer TPG Controller.

**class PfeifferTPGSerialCommunication(configuration)**

Bases: [hvl\\_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for Pfeiffer TPG controllers. Already predefines device-specific protocol parameters in config.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**query(cmd: str)** → str

Send a query, then read and returns the first line from the com port.

**Parameters cmd** – query message to send to the device

**Returns** first line read on the com

**Raises**

- [\*SerialCommunicationIOError\*](#) – when communication port is not opened
- [\*PfeifferTPGError\*](#) – if the device does not acknowledge the command or if the answer from the device is empty

**send\_command(cmd: str)** → None

Send a command to the device and check for acknowledgement.

**Parameters cmd** – command to send to the device

**Raises**

- [\*SerialCommunicationIOError\*](#) – when communication port is not opened
- [\*PfeifferTPGError\*](#) – if the answer from the device differs from the expected acknowledgement character ‘chr(6)’.

```
class PfeifferTPGSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str
= 'utf-8', encoding_error_handling: str =
'replace', wait_sec_read_text_nonempty:
Union[int, float] = 0.5, default_n_attempts_read_text_nonempty:
int = 10, port: Union[str, NoneType] =
None, baudrate: int = 9600, parity: Union[str,
hvl_ccb.comm.serial.SerialCommunicationParity]
= <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
hvl_ccb.comm.serial.SerialCommunicationStopbits]
= <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
hvl_ccb.comm.serial.SerialCommunicationBytesize]
= <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int,
float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

**baudrate:** `int = 9600`

Baudrate for Pfeiffer TPG controllers is 9600 baud

**bytesize:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8`  
One byte is eight bits long**force\_value** (`fieldname, value`)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**parity:** `Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'`

Pfeiffer TPG controllers do not use parity

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


**Returns** a list of strings containing all required keys.

**stopbits:** `Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1`  
Pfeiffer TPG controllers use one stop bit**terminator:** `bytes = b'\r\n'`

The terminator is <CR><LF>

```
timeout: Union[int, float] = 3
use 3 seconds timeout as default
```

## [hvl\\_ccb.dev.rs\\_rto1024](#)

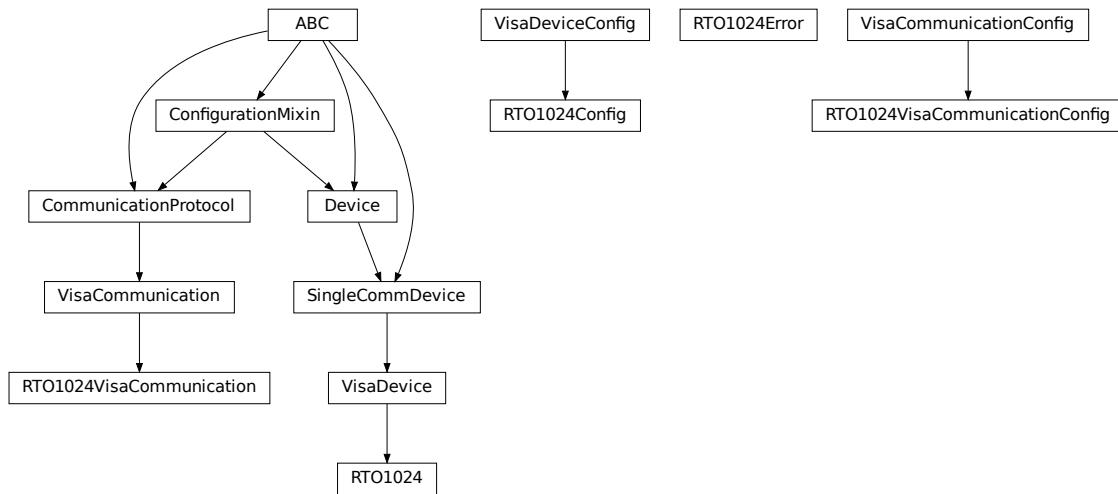

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

```
class RTO1024 (com: Union[hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunication,
                           hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunicationConfig, dict], dev_config:
                           Union[hvl_ccb.dev.rs_rto1024.RTO1024Config, dict])
Bases: hvl_ccb.dev.visa.VisaDevice
```

Device class for the Rhode & Schwarz RTO 1024 oscilloscope.

```
class TriggerModes (value=<object object>, names=None, module=None, type=None, start=1,
                     boundary=None)
Bases: hvl_ccb.utils.enum.AutoNumberNameEnum
```

Enumeration for the three available trigger modes.

**AUTO** = 1

**FREERUN** = 3

**NORMAL** = 2

**classmethod names()**

Returns a list of the available trigger modes. :return: list of strings

```
activate_measurements (meas_n: int, source: str, measurements: List[str], category: str = 'AMP-Time')
```

Activate the list of ‘measurements’ of the waveform ‘source’ in the measurement box number ‘meas\_n’. The list ‘measurements’ starts with the main measurement and continues with additional measurements of the same ‘category’.

### Parameters

- **meas\_n** – measurement number 1..8

- **source** – measurement source, for example C1W1
- **measurements** – list of measurements, the first one will be the main measurement.
- **category** – the category of measurements, by default AMPTime

**backup\_waveform** (*filename*: str) → None

Backup a waveform file from the standard directory specified in the device configuration to the standard backup destination specified in the device configuration. The filename has to be specified without .bin or path.

**Parameters** **filename** – The waveform filename without extension and path

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns** the VisaCommunication class

**file\_copy** (*source*: str, *destination*: str) → None

Copy a file from one destination to another on the oscilloscope drive. If the destination file already exists, it is overwritten without notice.

**Parameters**

- **source** – absolute path to the source file on the DSO filesystem
- **destination** – absolute path to the destination file on the DSO filesystem

**Raises** **RTO1024Error** – if the operation did not complete

**get\_acquire\_length()** → float

Gets the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 … 500 [s]
- Increment: 1E-12 [s]

**Returns** the time for one acquisition. Range: 250e-12 … 500 [s]

**get\_channel\_offset** (*channel*: int) → float

Gets the voltage offset of the indicated channel.

**Parameters** **channel** – is the channel number (1..4)

**Returns** channel offset voltage in V (value between -1 and 1)

**get\_channel\_position** (*channel*: int) → float

Gets the vertical position of the indicated channel.

**Parameters** **channel** – is the channel number (1..4)

**Returns** channel position in div (value between -5 and 5)

**get\_channel\_range** (*channel*: int) → float

Queries the channel range in V.

**Parameters** **channel** – is the input channel (1..4)

**Returns** channel range in V

**get\_channel\_scale** (*channel: int*) → float

Queries the channel scale in V/div.

**Parameters** **channel** – is the input channel (1..4)

**Returns** channel scale in V/div

**get\_channel\_state** (*channel: int*) → bool

Queries if the channel is active or not.

**Parameters** **channel** – is the input channel (1..4)

**Returns** True if active, else False

**get\_reference\_point** () → int

Gets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]

**Returns** the reference in %

**get\_repetitions** () → int

Get the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- \*RST = 1

**Returns** the number of waveforms to acquire

**get\_timestamps** () → List[float]

Gets the timestamps of all recorded frames in the history and returns them as a list of floats.

**Returns** list of timestamps in [s]

**Raises** *RTO1024Error* – if the timestamps are invalid

**list\_directory** (*path: str*) → List[Tuple[str, str, int]]

List the contents of a given directory on the oscilloscope filesystem.

**Parameters** **path** – is the path to a folder

**Returns** a list of filenames in the given folder

**load\_configuration** (*filename: str*) → None

Load current settings from a configuration file. The filename has to be specified without base directory and ‘.dfl’ extension.

**Information from the manual** *ReCaLL* calls up the instrument settings from an intermediate memory identified by the specified number. The instrument settings can be stored to this memory using the command \*SAV with the associated number. It also activates the instrument settings which are stored in a file and loaded using *MMEMory:LOAD:STATE*.

**Parameters** **filename** – is the name of the settings file without path and extension

**local\_display** (*state: bool*) → None

Enable or disable local display of the scope.

**Parameters** `state` – is the desired local display state

`prepare_ultra_segmentation() → None`

Make ready for a new acquisition in ultra segmentation mode. This function does one acquisition without ultra segmentation to clear the history and prepare for a new measurement.

`read_measurement(meas_n: int, name: str) → float`

**Parameters**

- `meas_n` – measurement number 1..8
- `name` – measurement name, for example “MAX”

**Returns** measured value

`run_continuous_acquisition() → None`

Start acquiring continuously.

`run_single_acquisition() → None`

Start a single or Nx acquisition.

`save_configuration(filename: str) → None`

Save the current oscilloscope settings to a file. The filename has to be specified without path and ‘.df1’ extension, the file will be saved to the configured settings directory.

**Information from the manual** `SAVe` stores the current instrument settings under the specified number in an intermediate memory. The settings can be recalled using the command `*RCL` with the associated number. To transfer the stored instrument settings to a file, use `MMEMory:STORe:STATe`.

**Parameters** `filename` – is the name of the settings file without path and extension

`save_waveform_history(filename: str, channel: int, waveform: int = 1) → None`

Save the history of one channel and one waveform to a .bin file. This function is used after an acquisition using sequence trigger mode (with or without ultra segmentation) was performed.

**Parameters**

- `filename` – is the name (without extension) of the file
- `channel` – is the channel number
- `waveform` – is the waveform number (typically 1)

**Raises** `RTO1024Error` – if storing waveform times out

`set_acquire_length(timerange: float) → None`

Defines the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 … 500 [s]
- Increment: 1E-12 [s]
- \*RST = 0.5 [s]

**Parameters** `timerange` – is the time for one acquisition. Range: 250e-12 … 500 [s]

`set_channel_offset(channel: int, offset: float) → None`

Sets the voltage offset of the indicated channel.

- Range: Dependent on the channel scale and coupling [V]
- Increment: Minimum 0.001 [V], may be higher depending on the channel scale and coupling
- \*RST = 0

### Parameters

- **channel1** – is the channel number (1..4)
- **offset** – Offset voltage. Positive values move the waveform down, negative values move it up.

**set\_channel\_position** (*channel: int, position: float*) → None

Sets the vertical position of the indicated channel as a graphical value.

- Range: -5.0 ... 5.0 [div]
- Increment: 0.02
- \*RST = 0

### Parameters

- **channel1** – is the channel number (1..4)
- **position** – is the position. Positive values move the waveform up, negative values move it down.

**set\_channel\_range** (*channel: int, v\_range: float*) → None

Sets the voltage range across the 10 vertical divisions of the diagram. Use the command alternatively instead of set\_channel\_scale.

- Range for range: Depends on attenuation factors and coupling. With 1:1 probe and external attenuations and 50 input coupling, the range is 10 mV to 10 V. For 1 M input coupling, it is 10 mV to 100 V. If the probe and/or external attenuation is changed, multiply the range values by the attenuation factors.
- Increment: 0.01
- \*RST = 0.5

### Parameters

- **channel1** – is the channel number (1..4)
- **v\_range** – is the vertical range [V]

**set\_channel\_scale** (*channel: int, scale: float*) → None

Sets the vertical scale for the indicated channel. The scale value is given in volts per division.

- Range for scale: depends on attenuation factor and coupling. With 1:1 probe and external attenuations and 50 input coupling, the vertical scale (input sensitivity) is 1 mV/div to 1 V/div. For 1 M input coupling, it is 1 mV/div to 10 V/div. If the probe and/or external attenuation is changed, multiply the values by the attenuation factors to get the actual scale range.
- Increment: 1e-3
- \*RST = 0.05

See also: set\_channel\_range

### Parameters

- **channel1** – is the channel number (1..4)
- **scale** – is the vertical scaling [V/div]

**set\_channel\_state** (*channel: int, state: bool*) → None

Switches the channel signal on or off.

**Parameters**

- **channel1** – is the input channel (1..4)
- **state** – is True for on, False for off

**set\_reference\_point** (*percentage: int*) → None

Sets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]
- \*RST = 50 [%]

**Parameters** **percentage** – is the reference in %

**set\_repetitions** (*number: int*) → None

Set the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- \*RST = 1

**Parameters** **number** – is the number of waveforms to acquire

**set\_trigger\_level** (*channel: int, level: float, event\_type: int = 1*) → None

Sets the trigger level for the specified event and source.

- Range: -10 to 10 V
- Increment: 1e-3 V
- \*RST = 0 V

**Parameters**

- **channel** – indicates the trigger source.
  - 1..4 = channel 1 to 4, available for all event types 1..3
  - 5 = external trigger input on the rear panel for analog signals, available for A-event type = 1
  - 6..9 = not available
- **level** – is the voltage for the trigger level in [V].
- **event\_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

**set\_trigger\_mode** (*mode: Union[str, hvl\_ccb.dev.rs\_rto1024.RTO1024.TriggerModes]*) → None

Sets the trigger mode which determines the behavior of the instrument if no trigger occurs.

**Parameters** **mode** – is either auto, normal, or freerun.

**Raises** **RTO1024Error** – if an invalid triggermode is selected

**set\_trigger\_source** (*channel: int, event\_type: int = 1*) → None

Set the trigger (Event A) source channel.

**Parameters**

- **channel1** – is the channel number (1..4)
- **event\_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

**start ()** → None

Start the RTO1024 oscilloscope and bring it into a defined state and remote mode.

**stop ()** → None

Stop the RTO1024 oscilloscope, reset events and close communication. Brings back the device to a state where local operation is possible.

**stop\_acquisition ()** → None

Stop any acquisition.

```
class RTO1024Config(waveforms_path: str, settings_path: str, backup_path: str, spoll_interval:
                      Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2, com-
                      mand_timeout_seconds: Union[int, float] = 60, wait_sec_short_pause:
                      Union[int, float] = 0.1, wait_sec_enable_history: Union[int, float] = 1,
                      wait_sec_post_acquisition_start: Union[int, float] = 2)
Bases: hvl_ccb.dev.visa.VisaDeviceConfig, hvl_ccb.dev.rs_rto1024._RTO1024ConfigDefaultsBase, hvl_ccb.dev.rs_rto1024._RTO1024ConfigBase
```

Configdataclass for the RTO1024 device.

**force\_value (fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys ()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults ()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys ()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**exception RTO1024Error**

Bases: Exception

**class RTO1024VisaCommunication (configuration)**

Bases: [hvl\\_ccb.comm.visa.VisaCommunication](#)

Specialization of VisaCommunication for the RTO1024 oscilloscope

**static config\_cls ()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class RTO1024VisaCommunicationConfig(host: str, interface_type: Union[str,
    hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]
    = <InterfaceType.TCPIP_INSTR: 2>, board: int = 0,
    port: int = 5025, timeout: int = 5000, chunk_size: int =
    204800, open_timeout: int = 1000, write_termination:
    str = '\n', read_termination: str = '\n', visa_backend:
    str = '')
```

Bases: `hvl_ccb.comm.visa.VisaCommunicationConfig`

Configuration dataclass for VisaCommunication with specifications for the RTO1024 device class.

**force\_value** (*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**interface\_type:** `Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]` =

Interface type of the VISA connection, being one of `InterfaceType`.

**classmethod keys()** → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → `Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.


**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → `Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

## `hvl_ccb.dev.se_il2t`

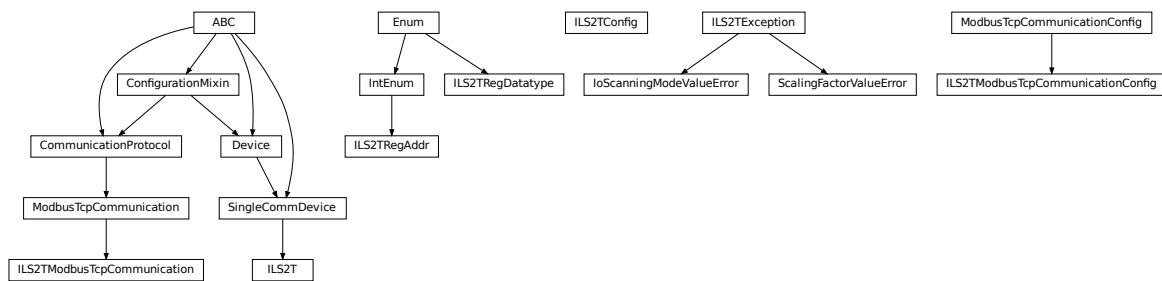

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

**class ILS2T** (*com*, *dev\_config=None*)

Bases: `hvl_ccb.dev.base.SingleCommDevice`

Schneider Electric ILS2T stepper drive class.

**ACTION\_JOG\_VALUE** = 0

The single action value for *ILS2T.Mode.JOG*

**class ActionsPtp**(*value*)

Bases: enum.IntEnum

Allowed actions in the point to point mode (*ILS2T.Mode.PTP*).

**ABSOLUTE\_POSITION** = 0

**RELATIVE\_POSITION\_MOTOR** = 2

**RELATIVE\_POSITION\_TARGET** = 1

**DEFAULT\_IO\_SCANNING\_CONTROL\_VALUES** = {'action': 2, 'continue\_after\_stop\_cu': 0, 'dis...

Default IO Scanning control mode values

**class Mode**(*value*)

Bases: enum.IntEnum

ILS2T device modes

**JOG** = 1

**PTP** = 3

**class Ref16Jog**(*value*)

Bases: enum.Flag

Allowed values for ILS2T ref\_16 register (the shown values are the integer representation of the bits), all in Jog mode = 1

**FAST** = 4

**NEG** = 2

**NEG\_FAST** = 6

**NONE** = 0

**POS** = 1

**POS\_FAST** = 5

**RegAddr**

Modbus Register Adresses

alias of [\*hvl\\_ccb.dev.se\\_ilis2t.ILS2TRegAddr\*](#)

**RegDatatype**

Modbus Register Datatypes

alias of [\*hvl\\_ccb.dev.se\\_ilis2t.ILS2TRegDatatype\*](#)

**class State**(*value*)

Bases: enum.IntEnum

State machine status values

**ON** = 6

**QUICKSTOP** = 7

**READY** = 4

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls()**

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**disable(log\_warn: bool = True, wait\_sec\_max: Optional[int] = None) → bool**

Disable the driver of the stepper motor and enable the brake.

Note: the driver cannot be disabled if the motor is still running.

**Parameters**

- **log\_warn** – if log a warning in case the motor cannot be disabled.
- **wait\_sec\_max** – maximal wait time for the motor to stop running and to disable it; by default, with *None*, use a config value

**Returns** *True* if disable request could and was sent, *False* otherwise.

**do\_ioscanning\_write(\*\*kwargs: int) → None**

Perform a write operation using IO Scanning mode.

**Parameters** **kwargs** – Keyword-argument list with options to send, remaining are taken from the defaults.

**enable() → None**

Enable the driver of the stepper motor and disable the brake.

**execute\_absolute\_position(position: int) → bool**

Execute a absolute position change, i.e. enable motor, perform absolute position change, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

**Parameters** **position** – absolute position of motor in user defined steps.

**Returns** *True* if actual position is as expected, *False* otherwise.

**execute\_relative\_step(steps: int) → bool**

Execute a relative step, i.e. enable motor, perform relative steps, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

**Parameters** **steps** – Number of steps.

**Returns** *True* if actual position is as expected, *False* otherwise.

**get\_dc\_volt() → float**

Read the DC supply voltage of the motor.

**Returns** DC input voltage.

**get\_error\_code() → Dict[int, Dict[str, Any]]**

Read all messages in fault memory. Will read the full error message and return the decoded values. At the end the fault memory of the motor will be deleted. In addition, reset\_error is called to re-enable the motor for operation.

**Returns** Dictionary with all information

**get\_position()** → int

Read the position of the drive and store into status.

**Returns** Position step value

**get\_status()** → Dict[str, int]

Perform an IO Scanning read and return the status of the motor.

**Returns** dict with status information.

**get\_temperature()** → int

Read the temperature of the motor.

**Returns** Temperature in degrees Celsius.

**jog\_run(direction: bool = True, fast: bool = False)** → None

Slowly turn the motor in positive direction.

**jog\_stop()** → None

Stop turning the motor in Jog mode.

**quickstop()** → None

Stops the motor with high deceleration rate and falls into error state. Reset with *reset\_error* to recover into normal state.

**reset\_error()** → None

Resets the motor into normal state after quick stop or another error occurred.

**set\_jog\_speed(slow: int = 60, fast: int = 180)** → None

Set the speed for jog mode. Default values correspond to startup values of the motor.

**Parameters**

- **slow** – RPM for slow jog mode.
- **fast** – RPM for fast jog mode.

**set\_max\_acceleration(rpm\_minute: int)** → None

Set the maximum acceleration of the motor.

**Parameters** **rpm\_minute** – revolution per minute per minute

**set\_max\_deceleration(rpm\_minute: int)** → None

Set the maximum deceleration of the motor.

**Parameters** **rpm\_minute** – revolution per minute per minute

**set\_max\_rpm(rpm: int)** → None

Set the maximum RPM.

**Parameters** **rpm** – revolution per minute ( 0 < rpm <= RPM\_MAX)

**Raises** *ILS2TException* – if RPM is out of range

**set\_ramp\_type(ramp\_type: int = -1)** → None

**Set the ramp type. There are two options available:** 0: linear ramp -1: motor optimized ramp

**Parameters** **ramp\_type** – 0: linear ramp | -1: motor optimized ramp

**start()** → None

Start this device.

**stop()** → None

Stop this device. Disables the motor (applies brake), disables access and closes the communication protocol.

**user\_steps** (*steps: int = 16384, revolutions: int = 1*) → None

Define steps per revolution. Default is 16384 steps per revolution. Maximum precision is 32768 steps per revolution.

#### Parameters

- **steps** – number of steps in *revolutions*.
- **revolutions** – number of revolutions corresponding to *steps*.

**write\_absolute\_position** (*position: int*) → None

Write instruction to turn the motor until it reaches the absolute position. This function does not enable or disable the motor automatically.

**Parameters** **position** – absolute position of motor in user defined steps.

**write\_relative\_step** (*steps: int*) → None

Write instruction to turn the motor the relative amount of steps. This function does not enable or disable the motor automatically.

**Parameters** **steps** – Number of steps to turn the motor.

```
class ILS2TConfig(rpm_max_init: numbers.Integral = 1500, wait_sec_post_enable: Union[int, float] = 1, wait_sec_max_disable: Union[int, float] = 10, wait_sec_post_cannot_disable: Union[int, float] = 1, wait_sec_post_relative_step: Union[int, float] = 2, wait_sec_post_absolute_position: Union[int, float] = 2)
```

Bases: object

Configuration for the ILS2T stepper motor device.

**clean\_values()****force\_value** (*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True****classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

```
rpm_max_init: numbers.Integral = 1500
    initial maximum RPM for the motor, can be set up to 3000 RPM. The user is allowed to set a new max
    RPM at runtime using ILS2T.set_max_rpm(), but the value must never exceed this configuration
    setting.

wait_sec_max_disable: Union[int, float] = 10
wait_sec_post_absolute_position: Union[int, float] = 2
wait_sec_post_CANNOT_DISABLE: Union[int, float] = 1
wait_sec_post_ENABLE: Union[int, float] = 1
wait_sec_post_relative_step: Union[int, float] = 2

exception ILS2TException
    Bases: Exception

    Exception to indicate problems with the SE ILS2T stepper motor.

class ILS2TModbusTcpCommunication(configuration)
    Bases: hvl_ccb.comm.modbus_tcp.ModbusTcpCommunication

    Specific implementation of Modbus/TCP for the Schneider Electric ILS2T stepper motor.

    static config_cls()
        Return the default configdataclass class.

        Returns a reference to the default configdataclass class

class ILS2TModbusTcpCommunicationConfig(host: str, unit: int = 255, port: int = 502)
    Bases: hvl_ccb.comm.modbus_tcp.ModbusTcpCommunicationConfig

    Configuration dataclass for Modbus/TCP communication specific for the Schneider Electric ILS2T stepper mo-
    tor.

    force_value(fieldname, value)
        Forces a value to a dataclass field despite the class being frozen.

        NOTE: you can define post_force_value method with same signature as this method to do extra processing
        after value has been forced on fieldname.

    Parameters
        • fieldname – name of the field
        • value – value to assign

    classmethod keys() → Sequence[str]
        Returns a list of all configdataclass fields key-names.

        Returns a list of strings containing all keys.

    classmethod optional_defaults() → Dict[str, object]
        Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
        on instantiation.

        Returns a list of strings containing all optional keys.

    classmethod required_keys() → Sequence[str]
        Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
        instantiation.

        Returns a list of strings containing all required keys.
```

```
unit: int = 255
The unit has to be 255 such that IO scanning mode works.
```

```
class ILS2TRegAddr (value)
```

Bases: enum.IntEnum

Modbus Register Adresses for Schneider Electric ILS2T stepper drive.

```
ACCESS_ENABLE = 282
```

```
FLT_INFO = 15362
```

```
FLT_MEM_DEL = 15112
```

```
FLT_MEM_RESET = 15114
```

```
IO_SCANNING = 6922
```

```
JOGN_FAST = 10506
```

```
JOGN_SLOW = 10504
```

```
POSITION = 7706
```

```
RAMP_ACC = 1556
```

```
RAMP_DECEL = 1558
```

```
RAMP_N_MAX = 1554
```

```
RAMP_TYPE = 1574
```

```
SCALE = 1550
```

```
TEMP = 7200
```

```
VOLT = 7198
```

```
class ILS2TRegDatatype (value=<object object>, names=None, module=None, type=None, start=1,
                           boundary=None)
```

Bases: aenum.Enum

Modbus Register Datatypes for Schneider Electric ILS2T stepper drive.

From the manual of the drive:

datatype	byte	min	max
INT8	1 Byte	-128	127
UINT8	1 Byte	0	255
INT16	2 Byte	-32_768	32_767
UINT16	2 Byte	0	65_535
INT32	4 Byte	-2_147_483_648	2_147_483_647
UINT32	4 Byte	0	4_294_967_295
BITS	just 32bits	N/A	N/A

```
INT32 = (-2147483648, 2147483647)
```

```
is_in_range (value: int) → bool
```

```
exception IoScanningModeValueError
```


Bases: *hvl\_ccb.dev.se\_ilst2t.ILS2TException*

Exception to indicate that the selected IO scanning mode is invalid.

```
exception ScalingFactorValueError
Bases: hvl\_ccb.dev.se\_ils2t.ILS2TException

Exception to indicate that a scaling factor value is invalid.
```

## [hvl\\_ccb.dev.sst\\_luminox](#)

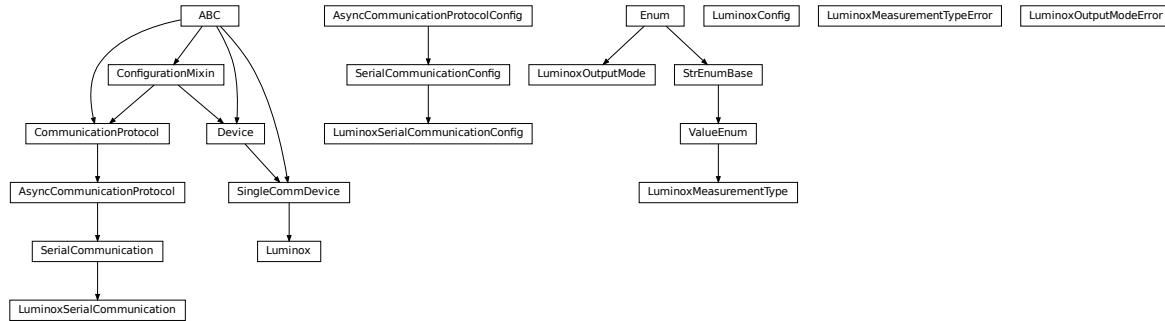

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.

Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

```
class Luminox(com, dev_config=None)
Bases: hvl\_ccb.dev.base.SingleCommDevice

Luminox oxygen sensor device class.

activate_output(mode: hvl_ccb.dev.sst_luminox.LuminoxOutputMode) → None
activate the selected output mode of the Luminox Sensor. :param mode: polling or streaming

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

query_polling(measurement: Union[str, hvl_ccb.dev.sst_luminox.LuminoxMeasurementType])
→ Union[Dict[Union[str, hvl_ccb.dev.sst_luminox.LuminoxMeasurementType], Union[float, int, str]], float, int, str]
Query a value or values of Luminox measurements in the polling mode, according to a given measurement type.

Parameters measurement – type of measurement

Returns value of requested measurement

Raises

- ValueError – when a wrong key for LuminoxMeasurementType is provided

```

- *LuminoxOutputModeError* – when polling mode is not activated
- *LuminoxMeasurementTypeError* – when expected measurement value is not read

**read\_streaming()** → Dict[Union[str, *hvl\_ccb.dev.sst\_luminox.LuminoxMeasurementType*], Union[float, int, str]]

Read values of Luminox in the streaming mode. Convert the single string into separate values.

**Returns** dictionary with *LuminoxMeasurementType.all\_measurements\_types()* keys and accordingly type-parsed values.

#### Raises

- *LuminoxOutputModeError* – when streaming mode is not activated
- *LuminoxMeasurementTypeError* – when any of expected measurement values is not read

**start()** → None

Start this device. Opens the communication protocol.

**stop()** → None

Stop the device. Closes also the communication protocol.

**class LuminoxConfig**(*wait\_sec\_post\_activate*: Union[int, float] = 0.5, *wait\_sec\_trials\_activate*: Union[int, float] = 0.1, *nr\_trials\_activate*: int = 5)

Bases: object

Configuration for the SST Luminox oxygen sensor.

**clean\_values()**

**force\_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass** = True

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**nr\_trials\_activate**: int = 5

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**wait\_sec\_post\_activate**: Union[int, float] = 0.5

```
    wait_sec_trials_activate: Union[int, float] = 0.1

class LuminoxMeasurementType(value=<object object>, names=None, module=None, type=None,
                                start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum
```

Measurement types for *LuminoxOutputMode.polling*.

The *all\_measurements* type will read values for the actual measurement types as given in *LuminoxOutputMode.all\_measurements\_types()*; it parses multiple single values using regexp's for other measurement types, therefore, no regexp is defined for this measurement type.

```
all_measurements = 'A'

classmethod all_measurements_types() → Tuple[hvl\_ccb.dev.sst\_luminox.LuminoxMeasurementType,
                                              ...]
A tuple of LuminoxMeasurementType enum instances which are actual measurements, i.e. not date of manufacture or software revision.

barometric_pressure = 'P'

property command

date_of_manufacture = '# 0'

parse_read_measurement_value(read_txt: str) → Union[Dict[Union[str,
                                                               hvl\_ccb.dev.sst\_luminox.LuminoxMeasurementType],
                                                               Union[float, int, str]], float, int, str]

partial_pressure_o2 = 'O'

percent_o2 = '%'

sensor_status = 'e'

serial_number = '# 1'

software_revision = '# 2'

temperature_sensor = 'T'
```

#### **LuminoxMeasurementTypeDict**

A typing hint for a dictionary holding *LuminoxMeasurementType* values. Keys are allowed as strings because *LuminoxMeasurementType* is of a *StrEnumBase* type.

alias of *Dict[Union[str, LuminoxMeasurementType], Union[float, int, str]]*

#### **exception LuminoxMeasurementTypeError**

Bases: *Exception*

Wrong measurement type for requested data

#### **LuminoxMeasurementTypeValue**

A typing hint for all possible *LuminoxMeasurementType* values as read in either streaming mode or in a polling mode with *LuminoxMeasurementType.all\_measurements*.

Beware: has to be manually kept in sync with *LuminoxMeasurementType* instances *cast\_type* attribute values.

alias of *Union[float, int, str]*

#### **class LuminoxOutputMode(value)**

Bases: *enum.Enum*

output mode.

**polling** = 1

**streaming** = 0

---

```
exception LuminoxOutputModeError
```

Bases: Exception

Wrong output mode for requested data

```
class LuminoxSerialCommunication(configuration)
```

Bases: `hvl_ccb.comm.serial.SerialCommunication`

Specific communication protocol implementation for the SST Luminox oxygen sensor. Already predefines device-specific protocol parameters in config.

```
static config_cls()
```

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

```
class LuminoxSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding_error_handling: str = 'replace', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

```
baudrate: int = 9600
```

Baudrate for SST Luminox is 9600 baud

```
bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
```

One byte is eight bits long

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

#### Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
classmethod keys() → Sequence[str]
```

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

```
classmethod optional_defaults() → Dict[str, object]
```

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

```

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'
    SST Luminox does not use parity

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.


stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
    SST Luminox does use one stop bit

terminator: bytes = b'\r\n'
    The terminator is CR LF

timeout: Union[int, float] = 3
    use 3 seconds timeout as default

```

## **hvl\_ccb.dev.technix**

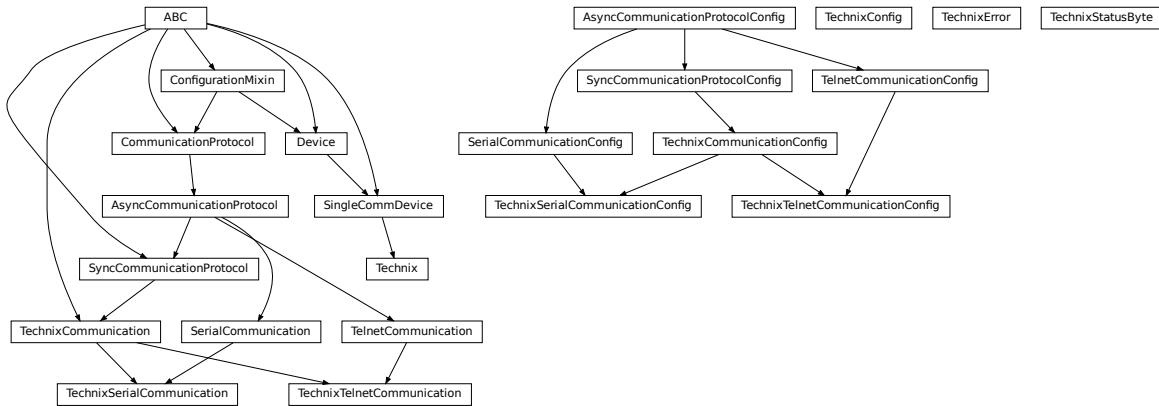

Device classes for “RS 232” and “Ethernet” Interfaces which are used to control power supplies from Technix. Manufacturer homepage: <https://www.technix-hv.com>

The Regulated power Supplies Series and Capacitor Chargers Series from Technix are series of low and high voltage direct current power supplies as well as capacitor chargers. The class *Technix* is tested with a CCR10KV-7,5KJ via an ethernet connection as well as a CCR15-P-2500-OP via a serial connection. Check the code carefully before using it with other devices or device series

**This Python package may support the following interfaces from Technix:**

- Remote Interface RS232
- Ethernet Remote Interface
- Optic Fiber Remote Interface

```

class Technix(com, dev_config)
    Bases: hvl_ccb.dev.base.SingleCommDevice

    static config_cls()
        Return the default configdataclass class.

```

**Returns** a reference to the default configdataclass class

```

property current
default_com_cls() → Union[Type[hvl_ccb.dev.technix.TechnixSerialCommunication],
                                Type[hvl_ccb.dev.technix.TechnixTelnetCommunication]]
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

get_status_byte() → hvl_ccb.dev.technix.TechnixStatusByte

property hv
property inhibit
property max_current
property max_voltage
property remote

start()
    Open the associated communication protocol.

stop()
    Close the associated communication protocol.

property voltage
property voltage_regulation

class TechnixCommunication(config)
    Bases: hvl_ccb.comm.base.SyncCommunicationProtocol, abc.ABC

    Generic communication class for Technix, which can be implemented via TechnixSerialCommunication or TechnixTelnetCommunication

query(command: str) → str
    Send a command to the interface and handle the status message. Eventually raises an exception.

        Parameters command – Command to send
        Raises TechnixError – if the connection is broken
        Returns Answer from the interface

class TechnixCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                    encoding_error_handling: str = 'replace',
                                    wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                    default_n_attempts_read_text_nonempty: int = 10)
    Bases: hvl_ccb.comm.base.SyncCommunicationProtocolConfig

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.

        Parameters
            • fieldname – name of the field
            • value – value to assign

classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

```

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**terminator: bytes = b'\r'**

The terminator is CR

**class TechnixConfig(communication\_channel: Union[Type[hvl\_ccb.dev.technix.TechnixSerialCommunication], Type[hvl\_ccb.dev.technix.TechnixTelnetCommunication]], max\_voltage: Union[int, float], max\_current: Union[int, float], polling\_interval\_sec: Union[int, float] = 4, post\_stop\_pause\_sec: Union[int, float] = 1, register\_pulse\_time: Union[int, float] = 0.1)**

Bases: object

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**communication\_channel: Union[Type[hvl\_ccb.dev.technix.TechnixSerialCommunication], Type[...]]**  
communication channel between computer and Technix

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

### Parameters

- **fieldname** – name of the field

- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**max\_current: Union[int, float]**

Maximal Output current

**max\_voltage: Union[int, float]**

Maximal Output voltage

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**polling\_interval\_sec: Union[int, float] = 4**

Polling interval in s to maintain to watchdog of the device

**post\_stop\_pause\_sec: Union[int, float] = 1**

Time to wait after stopping the device

```

register_pulse_time: Union[int, float] = 0.1
    Time for pulsing a register

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

exception TechnixError
    Bases: Exception

    Technix related errors.

class TechnixSerialCommunication(configuration)
    Bases: hvl_ccb.dev.technix.TechnixCommunication, hvl_ccb.comm.serial.SerialCommunication

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

class TechnixSerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8', encoding_error_handling: str = 'replace', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, float, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 2)
    Bases: hvl_ccb.dev.technix.TechnixCommunicationConfig, hvl_ccb.comm.serial.SerialCommunicationConfig

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.

Parameters

- fieldname – name of the field
- value – value to assign

classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

```

**classmethod required\_keys()** → Sequence[str]  
 Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

**class TechnixStatusByte (value: int)**

Bases: object

**msb\_first (idx: int)** → Optional[bool]

Give the Bit at position idx with MSB first

**Parameters** **idx** – Position of Bit as 1...8

**Returns**

**class TechnixTelnetCommunication (configuration)**

Bases: [hvl\\_ccb.comm.telnet.TelnetCommunication](#), [hvl\\_ccb.dev.technix.TechinxCommunication](#)

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**class TechnixTelnetCommunicationConfig (terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding\_error\_handling: str = 'replace', wait\_sec\_read\_text\_nonempty: Union[int, float] = 0.5, default\_n\_attempts\_read\_text\_nonempty: int = 10, host: Union[str, NoneType] = None, port: int = 4660, timeout: Union[int, float] = 0.2)**

Bases: [hvl\\_ccb.comm.telnet.TelnetCommunicationConfig](#), [hvl\\_ccb.dev.technix.TechinxCommunicationConfig](#)

**force\_value (fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**port: int = 4660**


Port at which Technix is listening

**classmethod required\_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.

[hvl\\_ccb.dev.tiepie](http://hvl_ccb.dev.tiepie)

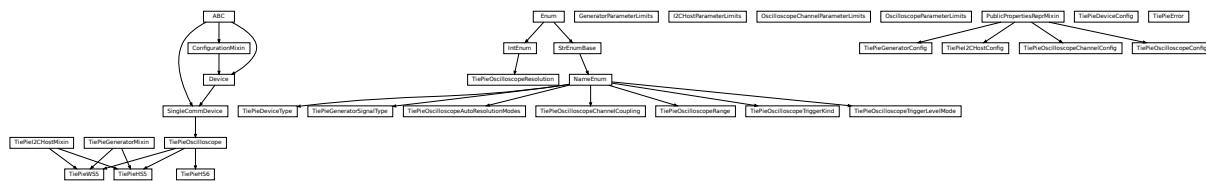

This module is a wrapper around libtiepie Oscilloscope devices; see <https://www.tiepie.com/en/libtiepie-sdk>.

The device classes adds simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and oscilloscope's channels. This includes extra validation and typing hints support.

To install libtiepie on Windows: The installation of the Python bindings “python-libtiepie” is done automatically with the dependencies of the hyl\_ccb. The additional DLL for Windows is included in that package.

On a Linux-system additional libraries have to be installed; see <https://www.tiepie.com/en/libtiepie-sdk/linux> .

On a Windows system, if you encounter an `OSError` like this:

```
    self._handle = _dlopen(self._name, mode)
OSError: [WinError 126] The specified module could not be found
```

most likely the python-libtiepie package was installed in your site-packages/ directory as a python-libtiepie-\*.egg file via python setup.py install or python setup.py develop command. In this case uninstall the library and re-install it using pip:

```
$ pip uninstall python-libtiepie  
$ pip install python-libtiepie
```

This should create `libtiepie/` folder. Alternatively, manually move the folder `libtiepie/` from inside of the `.egg` archive file to the containing it `site-packages/` directory (PyCharm's Project tool window supports reading and extracting from `.egg` archives).

```
class GeneratorParameterLimits (dev_gen: libtiepie.generator.Generator)
    Bases: object
```

#### Default limits for generator parameters

```
class I2CHostParameterLimits(dev_i2c: libtiepie.i2chost.I2CHost)
    Base: object
```

## Default limits for I2C host parameters

```
class OscilloscopeChannelParameterLimits(osc_channel: libtiepie.oscilloscopechannel.OscilloscopeChannel)
    Passes object
```

## Default limits for oscilloscope channel parameters

```
class OscilloscopeParameterLimits(dev_osc: libtiepie.oscilloscope.Oscilloscope):
    pass; object
```

## Default limits for oscilloscope parameters

```
class PublicPropertiesReprMixin
```

General purpose utility mixin that overwrites object representation to a one analogous to *dataclass* instances, but using public properties and their values instead of fields.

```
class TiePieDeviceConfig(serial_number: int, require_block_measurement_support: bool = True,
                         n_max_try_get_device: int = 10, wait_sec_retry_get_device: Union[int,
                         float] = 1.0, is_data_ready_polling_interval_sec: Union[int, float] =
                         0.01)
Bases: object
Configuration dataclass for TiePie

clean_values()
force_value(fieldname, value)
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define post_force_value method with same signature as this method to do extra processing
after value has been forced on fieldname.

Parameters

- fieldname – name of the field
- value – value to assign


is_configdataclass = True
is_data_ready_polling_interval_sec: Union[int, float] = 0.01
classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

n_max_try_get_device: int = 10
classmethod optional_defaults() → Dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
on instantiation.

Returns a list of strings containing all optional keys.

require_block_measurement_support: bool = True
classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
instantiation.

Returns a list of strings containing all required keys.

serial_number: int
wait_sec_retry_get_device: Union[int, float] = 1.0
class TiePieDeviceType(value=<object object>, names=None, module=None, type=None, start=1,
                      boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
TiePie device type.

GENERATOR = 2
I2C = 4
OSCILLOSCOPE = 1

exception TiePieError
Bases: Exception
Error of the class TiePie
```

```

class TiePieGeneratorConfig(dev_gen: libtiepie.generator.Generator)
Bases: hvl_ccb.dev.tiepie.PublicPropertiesReprMixin
Generator's configuration with cleaning of values in properties setters.

property amplitude
clean_amplitude(amplitude: float) → float
static clean_enabled(enabled: bool) → bool
clean_frequency(frequency: float) → float
clean_offset(offset: float) → float
static clean_signal_type(signal_type: Union[int, hvl_ccb.dev.tiepie.TiePieGeneratorSignalType])
→ hvl_ccb.dev.tiepie.TiePieGeneratorSignalType

property enabled
property frequency
property offset
property signal_type

class TiePieGeneratorMixin(com, dev_config)
Bases: object
TiePie Generator sub-device.

A wrapper for the libtiepie.generator.Generator class. To be mixed in with TiePieOscilloscope base class.

config_gen: Optional[hvl_ccb.dev.tiepie.TiePieGeneratorConfig]
Generator's dynamical configuration.

generator_start()
Start signal generation.

generator_stop()
Stop signal generation.

start() → None
Start the Generator.

stop() → None
Stop the generator.

class TiePieGeneratorSignalType(value=<object object>, names=None, module=None,
type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
An enumeration.

ARBITRARY = 32
DC = 8
NOISE = 16
PULSE = 64
SINE = 1
SQUARE = 4
TRIANGLE = 2
UNKNOWN = 0

```

```
class TiePieHS5(com, dev_config)
Bases: hvl_ccb.dev.tiepie.TiePieI2CHostMixin, hvl_ccb.dev.tiepie.TiePieGeneratorMixin, hvl_ccb.dev.tiepie.TiePieOscilloscope
TiePie HS5 device.

config_gen: Optional[hvl_ccb.dev.tiepie.TiePieGeneratorConfig]
Generator's dynamical configuration.

class TiePieHS6(com, dev_config)
Bases: hvl_ccb.dev.tiepie.TiePieOscilloscope
TiePie HS6 DIFF device.

class TiePieI2CHostConfig(dev_i2c: libtiepie.i2chost.I2CHost)
Bases: hvl_ccb.dev.tiepie.PublicPropertiesReprMixin
I2C Host's configuration with cleaning of values in properties setters.

class TiePieI2CHostMixin(com, dev_config)
Bases: object
TiePie I2CHost sub-device.

A wrapper for the libtiepie.i2chost.I2CHost class. To be mixed in with TiePieOscilloscope base class.

config_i2c: Optional[hvl_ccb.dev.tiepie.TiePieI2CHostConfig]
I2C host's dynamical configuration.

start() → None
Start the I2C Host.

stop() → None
Stop the I2C host.

class TiePieOscilloscope(com, dev_config)
Bases: hvl_ccb.dev.base.SingleCommDevice
TiePie oscilloscope.

A wrapper for TiePie oscilloscopes, based on the class libtiepie.oscilloscope.Oscilloscope with simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and its channels, including extra validation and typing hints support for configurations.

Note that, in contrast to libtiepie library, since all physical TiePie devices include an oscilloscope, this is the base class for all physical TiePie devices. The additional TiePie sub-devices: "Generator" and "I2CHost", are mixed-in to this base class in subclasses.

The channels use I..N numbering (not 0..N-1), as in, e.g., the Multi Channel software.

property channels_enabled
Yield numbers of enabled channels.

Returns Numbers of enabled channels

collect_measurement_data(timeout: Optional[Union[int, float]] = 0) → Optional[numumpy.ndarray]
Try to collect the data from TiePie; return None if data is not ready.

Parameters timeout – The timeout to wait until data is available. This option makes this function blocking the code. timeout = None blocks the code infinitely till data will be available. Per default, the timeout is set to 0: The function will not block.

Returns Measurement data of only enabled channels and time vector in a 2D-numpy.ndarray with float sample data; or None if there is no data available.
```

**static config\_cls()** → Type[hvl\_ccb.dev.tiepie.TiePieDeviceConfig]

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**config\_osc:** Optional[hvl\_ccb.dev.tiepie.TiePieOscilloscopeConfig]

Oscilloscope's dynamical configuration.

**config\_osc\_channel\_dict:** Dict[int, hvl\_ccb.dev.tiepie.TiePieOscilloscopeChannelConfig]

Channel configuration. A *dict* mapping actual channel number, numbered *I..N*, to channel configuration.

The channel info is dynamically read from the device only on the first *start()*; beforehand the *dict* is empty.

**static default\_com\_cls()** → Type[hvl\_ccb.comm.base.NullCommunicationProtocol]

Get the class for the default communication protocol used with this device.

**Returns** the type of the standard communication protocol for this device

**force\_trigger()** → None

Forces the TiePie to trigger with a software sided trigger event.

**Return None**

**Raises TiePieError** – when device is not started or status of underlying device gives an error

**is\_measurement\_data\_ready()** → bool

Reports if TiePie has data which is ready to collect

**Returns** if the data is ready to collect.

**Raises TiePieError** – when device is not started or status of underlying device gives an error

**is\_triggered()** → bool

Reports if TiePie has triggered. Maybe data is not yet available. One can check with the function *is\_measurement\_data\_ready()*.

**Returns** if a trigger event occurred

**static list\_devices()** → libtiepie.devicelist.DeviceList

List available TiePie devices.

**Returns** libtiepie up to date list of devices

**property n\_channels**

Number of channels in the oscilloscope.

**Returns** Number of channels.

**start()** → None

Start the oscilloscope.

**start\_measurement()** → None

Start a measurement using set configuration.

**Raises TiePieError** – when device is not started or status of underlying device gives an error

**stop()** → None

Stop the oscilloscope.

**class TiePieOscilloscopeAutoResolutionModes** (value=<object object>, names=None, module=None, type=None, start=1, boundary=None)

Bases: hvl\_ccb.utils.enum.NameEnum

An enumeration.

**ALL = 4**

```
DISABLED = 1
NATIVEONLY = 2
UNKNOWN = 0

class TiePieOscilloscopeChannelConfig(ch_number: int, channel: libtiepie.oscilloscopechannel.OscilloscopeChannel)
Bases: hvl_ccb.dev.tiepie.PublicPropertiesReprMixin

Oscilloscope's channel configuration, with cleaning of values in properties setters as well as setting and reading them on and from the device's channel.

static clean_coupling(coupling: Union[str, hvl_ccb.dev.tiepie.TiePieOscilloscopeChannelCoupling]) → hvl_ccb.dev.tiepie.TiePieOscilloscopeChannelCoupling
static clean_enabled(enabled: bool) → bool
clean_input_range(input_range: Union[float, hvl_ccb.dev.tiepie.TiePieOscilloscopeRange]) → hvl_ccb.dev.tiepie.TiePieOscilloscopeRange
clean_probe_offset(probe_offset: float) → float
static clean_safe_ground_enabled(safe_ground_enabled: bool) → bool
static clean_trigger_enabled(trigger_enabled)
clean_trigger_hysteresis(trigger_hysteresis: float) → float
static clean_trigger_kind(trigger_kind: Union[str, hvl_ccb.dev.tiepie.TiePieOscilloscopeTriggerKind]) → hvl_ccb.dev.tiepie.TiePieOscilloscopeTriggerKind
clean_trigger_level(trigger_level: Union[int, float]) → float
static clean_trigger_level_mode(level_mode: Union[str, hvl_ccb.dev.tiepie.TiePieOscilloscopeTriggerLevelMode]) → hvl_ccb.dev.tiepie.TiePieOscilloscopeTriggerLevelMode

property coupling
property enabled
property has_safe_ground
    Check whether bound oscilloscope device has "safe ground" option
        Returns bool: 1=safe ground available
property input_range
property probe_offset
property safe_ground_enabled
property trigger_enabled
property trigger_hysteresis
property trigger_kind
property trigger_level
property trigger_level_mode

class TiePieOscilloscopeChannelCoupling(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ACA = 8
```

```

ACV = 2
DCA = 4
DCV = 1

class TiePieOscilloscopeConfig(dev_osc: libtiepie.oscilloscope.Oscilloscope)
Bases: hvl_ccb.dev.tiepie.PublicPropertiesReprMixin

Oscilloscope's configuration with cleaning of values in properties setters.

property auto_resolution_mode
static clean_auto_resolution_mode(auto_resolution_mode: Union[int,  

                                         hvl_ccb.dev.tiepie.TiePieOscilloscopeAutoResolutionModes])  

                                         → hvl_ccb.dev.tiepie.TiePieOscilloscopeAutoResolutionModes

clean_pre_sample_ratio(pre_sample_ratio: float) → float
clean_record_length(record_length: Union[int, float]) → int
static clean_resolution(resolution: Union[int, hvl_ccb.dev.tiepie.TiePieOscilloscopeResolution])  

                                         → hvl_ccb.dev.tiepie.TiePieOscilloscopeResolution

clean_sample_frequency(sample_frequency: float) → float
clean_trigger_timeout(trigger_timeout: Optional[Union[int, float]]) → float

property pre_sample_ratio
property record_length
property resolution
property sample_frequency
property trigger_timeout

class TiePieOscilloscopeRange(value=<object object>, names=None, module=None, type=None,  

                                 start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

EIGHTY_VOLT = 80
EIGHT_HUNDRED_MILLI_VOLT = 0.8
EIGHT_VOLT = 8
FORTY_VOLT = 40
FOUR_HUNDRED_MILLI_VOLT = 0.4
FOUR_VOLT = 4
TWENTY_VOLT = 20
TWO_HUNDRED_MILLI_VOLT = 0.2
TWO_VOLT = 2

static suitable_range(value)

class TiePieOscilloscopeResolution(value=<object object>, names=None, module=None,  

                                         type=None, start=1, boundary=None)
Bases: aenum.IntEnum

An enumeration.

```

```
EIGHT_BIT = 8
FOURTEEN_BIT = 14
SIXTEEN_BIT = 16
TWELVE_BIT = 12

class TiePieOscilloscopeTriggerKind(value=<object object>, names=None, module=None,
                                      type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ANY = 16
FALLING = 2
RISING = 1
RISING_OR_FALLING = 16

class TiePieOscilloscopeTriggerLevelMode(value=<object object>, names=None, mod-
                                           ule=None, type=None, start=1, bound-
                                           ary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ABSOLUTE = 2
RELATIVE = 1
UNKNOWN = 0

class TiePieWS5(com, dev_config)
Bases:           hvl_ccb.dev.tiepie.TiePieI2CHostMixin,      hvl_ccb.dev.tiepie.
               TiePieGeneratorMixin, hvl_ccb.dev.tiepie.TiePieOscilloscope

TiePie WS5 device.

config_i2c: Optional[hvl_ccb.dev.tiepie.TiePieI2CHostConfig]
I2C host's dynamical configuration.

get_device_by_serial_number(serial_number: int, device_type: Union[str, Tuple[int, _LtpDeviceReturnType]], n_max_try_get_device: int = 10, wait_sec_retry_get_device: float = 1.0) → _LtpDeviceReturn-Type
Open and return handle of TiePie device with a given serial number
```

#### Parameters

- **serial\_number** – int serial number of the device
- **device\_type** – a *TiePieDeviceType* instance containing device identifier (int number) and its corresponding class, both from *libtiepie*, or a string name of such instance
- **n\_max\_try\_get\_device** – maximal number of device list updates (int number)
- **wait\_sec\_retry\_get\_device** – waiting time in seconds between retries (int number)

**Returns** Instance of a *libtiepie* device class according to the specified *device\_type*

#### Raises

- ***TiePieError*** – when there is no device with given serial number

- **ValueError** – when *device\_type* is not an instance of *TiePieDeviceType*

**log\_set** (*prop\_name*: str, *prop\_value*: object, *value\_suffix*: str = '') → None

**wrap\_libtiepie\_exception** (*func*: Callable) → Callable

Decorator wrapper for *libtiepie* methods that use *libtiepie.library.check\_last\_status\_raise\_on\_error()* calls.

**Parameters** **func** – Function or method to be wrapped

**Raises** **TiePieError** – instead of *LibTiePieException* or one of its subtypes.

**Returns** whatever *func* returns

## hvl\_ccb.dev.utils

### Poller

**class Poller** (*spoll\_handler*: Callable, *polling\_delay\_sec*: Union[int, float] = 0, *polling\_interval\_sec*: Union[int, float] = 1, *polling\_timeout\_sec*: Optional[Union[int, float]] = None)

Bases: object

Poller class wrapping *concurrent.futures.ThreadPoolExecutor* which enables passing of results and errors out of the polling thread.

**is\_polling** () → bool

Check if device status is being polled.

**Returns** *True* when polling thread is set and alive

**start\_polling** () → bool

Start polling.

**Returns** *True* if was not polling before, *False* otherwise

**stop\_polling** () → bool

Stop polling.

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

**Returns** *True* if was polling before, *False* otherwise, and last result of the polling function call.

**Raises** polling function exceptions


**wait\_for\_polling\_result** ()

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

**Returns** polling function result

**Raises** polling function errors

## hvl\_ccb.dev.visa

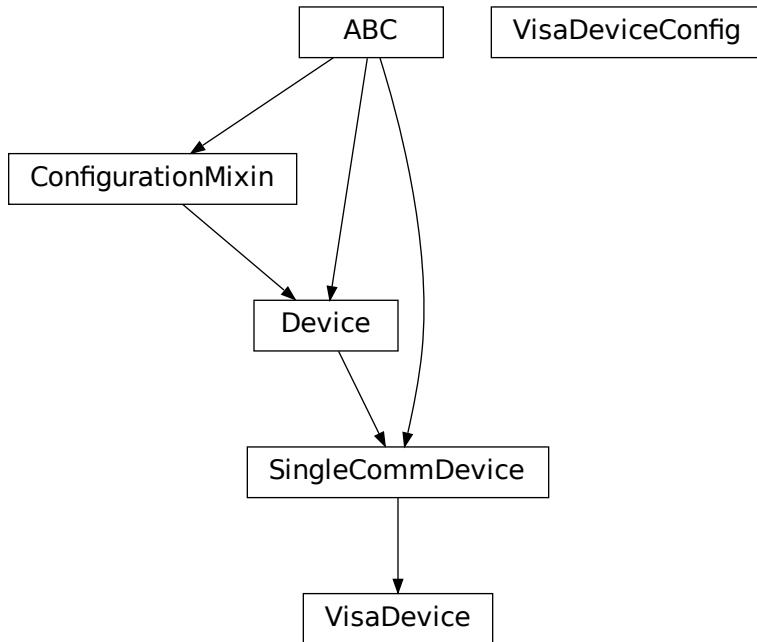

```
class VisaDevice(com: Union[hvl_ccb.comm.visa.VisaCommunication,
                           hvl_ccb.comm.visa.VisaCommunicationConfig, dict], dev_config: Optional[Union[hvl_ccb.dev.visa.VisaDeviceConfig, dict]] = None)
Bases: hvl_ccb.dev.base.SingleCommDevice
```

Device communicating over the VISA protocol using VisaCommunication.

**static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**static default\_com\_cls() → Type[hvl\_ccb.comm.visa.VisaCommunication]**

Return the default communication protocol for this device type, which is VisaCommunication.

**Returns** the VisaCommunication class

**get\_error\_queue() → str**

Read out error queue and logs the error.

**Returns** Error string

**get\_identification() → str**

Queries “\*IDN?” and returns the identification string of the connected device.

**Returns** the identification string of the connected device

**reset() → None**

Send “\*RST” and “\*CLS” to the device. Typically sets a defined state.

**spoll\_handler()**

Reads the status byte and decodes it. The status byte STB is defined in IEEE 488.2. It provides a rough overview of the instrument status.

**Returns****start() → None**

Start the VisaDevice. Sets up the status poller and starts it.

**Returns****stop() → None**

Stop the VisaDevice. Stops the polling thread and closes the communication protocol.

**Returns****wait\_operation\_complete(timeout: Optional[float] = None) → bool**

Waits for a operation complete event. Returns after timeout [s] has expired or the operation complete event has been caught.

**Parameters** **timeout** – Time in seconds to wait for the event; *None* for no timeout.

**Returns** True, if OPC event is caught, False if timeout expired

**class VisaDeviceConfig(spoll\_interval: Union[int, float] = 0.5, spoll\_start\_delay: Union[int, float] = 2)**  
Bases: hvl\_ccb.dev.visa.\_VisaDeviceConfigDefaultsBase, hvl\_ccb.dev.visa.\_VisaDeviceConfigBase

Configdataclass for a VISA device.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


## Module contents

Devices subpackage.

### [hvl\\_ccb.utils](#)

#### Submodules

##### [hvl\\_ccb.utils.enum](#)

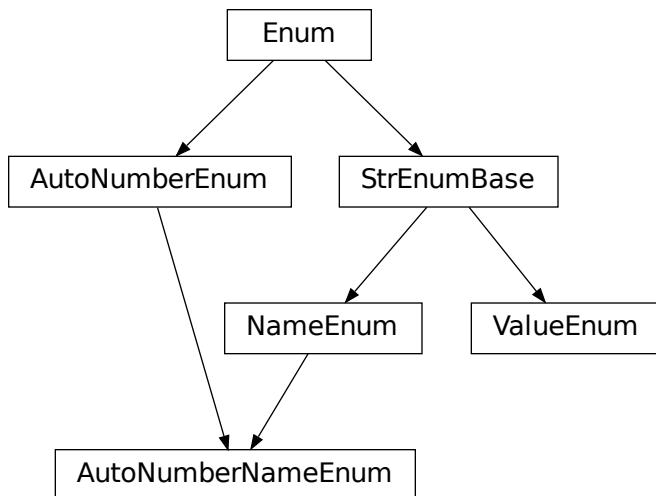

```
class AutoNumberNameEnum(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.NameEnum, aenum.AutoNumberEnum
```

Auto-numbered enum with names used as string representation, and with lookup and equality based on this representation.

```
class NameEnum(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.StrEnumBase
```

Enum with names used as string representation, and with lookup and equality based on this representation.

```
class StrEnumBase(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.Enum
```

String representation-based equality and lookup.

```
class ValueEnum(value=<object object>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.StrEnumBase
```

Enum with string representation of values used as string representation, and with lookup and equality based on this representation.

Attention: to avoid errors, best use together with *unique* enum decorator.

## **hvl\_ccb.utils.typing**

Additional Python typing module utilities

### **Number**

Typing hint auxiliary for a Python base number types: *int* or *float*.

alias of Union[int, float]

### **check\_generic\_type (value, type\_, name='instance')**

Check if *value* is of a generic type *type\_*. Raises *TypeError* if it's not.

#### **Parameters**

- **name** – name to report in case of an error
- **value** – value to check
- **type** – generic type to check against

### **is\_generic\_type\_hint (type\_)**


Check if class is a generic type, for example *Union[int, float]*, *List[int]*, or *List*.

#### **Parameters **type** – type to check**

## **Module contents**

### **4.1.2 Submodules**

## **hvl\_ccb.configuration**

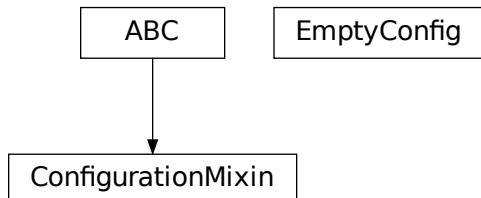

Facilities providing classes for handling configuration for communication protocols and devices.

**class ConfigurationMixin(configuration)**  
Bases: abc.ABC

Mixin providing configuration to a class.

**property config**  
ConfigDataclass property.

**Returns** the configuration

**abstract static config\_cls()**

Return the default configdataclass class.

**Returns** a reference to the default configdataclass class

**configuration\_save\_json(path: str) → None**

Save current configuration as JSON file.

**Parameters** **path** – path to the JSON file.

**classmethod from\_json(filename: str)**

Instantiate communication protocol using configuration from a JSON file.

**Parameters** **filename** – Path and filename to the JSON configuration

**class EmptyConfig**

Bases: object

Empty configuration dataclass.

**clean\_values()**

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

**force\_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post\_force\_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

**is\_configdataclass = True**

**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

**Returns** a list of strings containing all keys.

**classmethod optional\_defaults() → Dict[str, object]**

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

**Returns** a list of strings containing all optional keys.

**classmethod required\_keys() → Sequence[str]**

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

**Returns** a list of strings containing all required keys.


**configdataclass(direct\_decoration=None, frozen=True)**

Decorator to make a class a configdataclass. Types in these dataclasses are enforced. Implement a function *clean\_values(self)* to do additional checking on value ranges etc.

It is possible to inherit from a configdataclass and re-decorate it with @configdataclass. In a subclass, default values can be added to existing fields. Note: adding additional non-default fields is prone to errors, since the order has to be respected through the whole chain (first non-default fields, only then default-fields).

**Parameters** `frozen` – defaults to True. False allows to later change configuration values. Attention: if configdataclass is not frozen and a value is changed, typing is not enforced anymore!

## `hvl_ccb.experiment_manager`

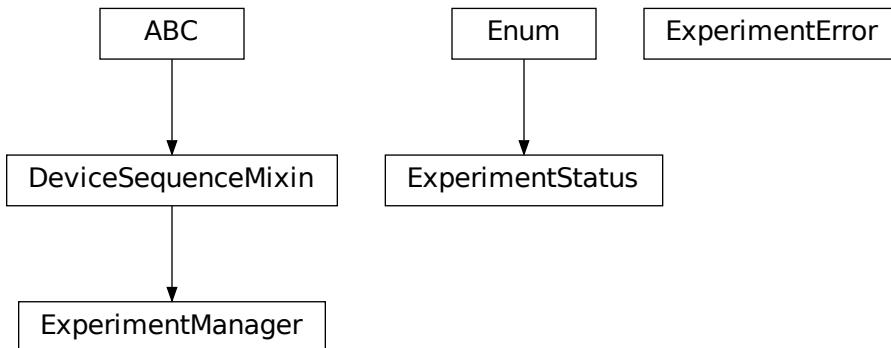

Main module containing the top level `ExperimentManager` class. Inherit from this class to implement your own experiment functionality in another project and it will help you start, stop and manage your devices.

### `exception ExperimentError`

Bases: `Exception`

Exception to indicate that the current status of the experiment manager is on ERROR and thus no operations can be made until reset.

### `class ExperimentManager(*args, **kwargs)`

Bases: `hvl_ccb.dev.base.DeviceSequenceMixin`

Experiment Manager can start and stop communication protocols and devices. It provides methods to queue commands to devices and collect results.

#### `add_device(name: str, device: hvl_ccb.dev.base.Device) → None`

Add a new device to the manager. If the experiment is running, automatically start the device. If a device with this name already exists, raise an exception.

##### Parameters

- `name` – is the name of the device.
- `device` – is the instantiated `Device` object.

**Raises** `DeviceExistingException` –

#### `devices_failed_start: Dict[str, hvl_ccb.dev.base.Device]`

Dictionary of named device instances from the sequence for which the most recent `start()` attempt failed.

Empty if `stop()` was called last; cf. `devices_failed_stop`.

#### `devices_failed_stop: Dict[str, hvl_ccb.dev.base.Device]`

Dictionary of named device instances from the sequence for which the most recent `stop()` attempt failed.

Empty if `start()` was called last; cf. `devices_failed_start`.

```
finish() → None
    Stop experimental setup, stop all devices.

is_error() → bool
    Returns true, if the status of the experiment manager is error.
    Returns True if on error, false otherwise

is_finished() → bool
    Returns true, if the status of the experiment manager is finished.
    Returns True if finished, false otherwise

is_running() → bool
    Returns true, if the status of the experiment manager is running.
    Returns True if running, false otherwise

run() → None
    Start experimental setup, start all devices.

start() → None
    Alias for ExperimentManager.run()

property status
    Get experiment status.
    Returns experiment status enum code.

stop() → None
    Alias for ExperimentManager.finish()

class ExperimentStatus (value)
    Bases: enum.Enum
    Enumeration for the experiment status
        ERROR = 5
        FINISHED = 4
        FINISHING = 3
        INITIALIZED = 0
        INITIALIZING = -1
        RUNNING = 2
        STARTING = 1
```

#### 4.1.3 Module contents

Top-level package for HVL Common Code Base.

## **CONTRIBUTING**

Contributions are welcome, and they are greatly appreciated! Every little bit helps, and credit will always be given.

You can contribute in many ways:

### **5.1 Types of Contributions**

#### **5.1.1 Report Bugs**

Report bugs at [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/issues](https://gitlab.com/ethz_hvl/hvl_ccb/issues).

If you are reporting a bug, please include:

- Your operating system name and version.
- Any details about your local setup that might be helpful in troubleshooting.
- Detailed steps to reproduce the bug.

#### **5.1.2 Fix Bugs**

Look through the GitLab issues for bugs. Anything tagged with “bug” and “help wanted” is open to whoever wants to implement it.

#### **5.1.3 Implement Features**

Look through the GitLab issues for features. Anything tagged with “enhancement” and “help wanted” is open to whoever wants to implement it.

#### **5.1.4 Write Documentation**

HVL Common Code Base could always use more documentation, whether as part of the official HVL Common Code Base docs, in docstrings, or even on the web in blog posts, articles, and such.

### 5.1.5 Submit Feedback

The best way to send feedback is to file an issue at [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/issues](https://gitlab.com/ethz_hvl/hvl_ccb/issues).

If you are proposing a feature:

- Explain in detail how it would work.
- Keep the scope as narrow as possible, to make it easier to implement.
- Remember that this is a volunteer-driven project, and that contributions are welcome :)

## 5.2 Get Started!

Ready to contribute? Here's how to set up *hvl\_ccb* for local development.

1. Clone *hvl\_ccb* repo from GitLab.

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

2. Install your local copy into a virtualenv. Assuming you have `virtualenvwrapper` installed, this is how you set up your fork for local development:

```
$ mkvirtualenv hvl_ccb
$ cd hvl_ccb/
$ pip install -e .[all]
$ pip install -r requirements_dev.txt
```

3. Create a branch for local development:

```
$ git checkout -b name-of-your-bugfix-or-feature
```

Now you can make your changes locally.

4. When you're done making changes, check that your changes pass flake8 and the tests, including testing other Python versions with tox:

```
$ flake8 hvl_ccb tests
$ python setup.py test or py.test
$ tox
```

To get flake8 and tox, just pip install them into your virtualenv. You can also use the provided make-like shell script to run flake8 and tests:

```
$ ./make.sh lint
$ ./make.sh test
```

5. Commit your changes and push your branch to GitLab:

```
$ git add .
$ git commit -m "Your detailed description of your changes."
$ git push origin name-of-your-bugfix-or-feature
```

6. Submit a merge request through the GitLab website.

## 5.3 Merge Request Guidelines

Before you submit a merge request, check that it meets these guidelines:

1. The merge request should include tests.
2. If the merge request adds functionality, the docs should be updated. Put your new functionality into a function with a docstring, and add the feature to the list in README.rst.
3. The merge request should work for Python 3.7. Check [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/merge\\_requests](https://gitlab.com/ethz_hvl/hvl_ccb/merge_requests) and make sure that the tests pass for all supported Python versions.

## 5.4 Tips

- To run tests from a single file:

```
$ py.test tests/test_hvl_ccb.py
```

or a single test function:

```
$ py.test tests/test_hvl_ccb.py::test_command_line_interface
```

- If your tests are slow, profile them using the pytest-profiling plugin:

```
$ py.test tests/test_hvl_ccb.py --profile
```

or for a graphical overview (you need a SVG image viewer):

```
$ py.test tests/test_hvl_ccb.py --profile-svg
$ open prof/combined.svg
```

- To add dependency, edit appropriate `*requirements` variable in the `setup.py` file and re-run:

```
$ python setup.py develop
```

- To generate a PDF version of the Sphinx documentation instead of HTML use:

```
$ rm -rf docs/hvl_ccb.rst docs/modules.rst docs/_build && sphinx-apidoc -o docs/
  ↪hvl_ccb && python -msphinx -M latexpdf docs/ docs/_build
```

This command can also be run through the make-like shell script:

```
$ ./make.sh docs-pdf
```

This requires a local installation of a LaTeX distribution, e.g. MikTeX.

## 5.5 Deploying

A reminder for the maintainers on how to deploy. Create `release-N.M.K` branch. Make sure all your changes are committed. Update or create entry in `HISTORY.rst` file, update features list in `README.rst` file and update API docs:

```
$ make docs
```

Commit all of the above and then run:

```
$ bumpversion patch # possible: major / minor / patch
$ git push
$ git push --tags
$ make release
```

Merge the release branch into master and devel branches with `--no-ff` flag and delete the release branch:

```
$ git checkout master
$ git merge --no-ff release-N.M.K
$ git checkout devel
$ git merge --no-ff release-N.M.K
$ git push --delete origin release-N.M.K
$ git branch --delete release-N.M.K
```

Finally, go to [https://gitlab.com/ethz\\_hvl/hvl\\_ccb/tags/](https://gitlab.com/ethz_hvl/hvl_ccb/tags/), select the latest release tag, press “Edit release notes” and add release notes (corresponding entry from `HISTORY.rst` file, but consider also additional brief header or synopsis if needed).

## CREDITS

### 6.1 Maintainers

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>

### 6.2 Authors

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- David Gruber <dev@davidgruber.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Alise Chachereau <chachereau@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>

### 6.3 Contributors

- Luca Nembrini <lucane@student.ethz.ch>
- Maria Pukhlyakova <mariia.pukhliakova@id.ethz.ch>
- Raphael Faerber <raphael.faerber@eeh.ee.ethz.ch>
- Ruben Stadler <rstadler@student.ethz.ch>



## HISTORY

### 7.1 0.6.1 (2021-05-08)

- In `dev.tiepie`: \* dynamically set oscilloscope's channel limits in `OscilloscopeChannelParameterLimits`: `input_range` and `trigger_level_abs`, incl. update of latter on each change of `input_range` value of a `TiePieOscilloscopeChannelConfig` instances;
  - quick fix for opening of combined instruments by disabling `OscilloscopeParameterLimits.trigger_delay` (an advanced feature);
  - enable automatic devices detection to be able to find network devices with `TiePieOscilloscope.list_devices()`.
- Fix `examples/example_labjack.py`.
- Improved logging: consistently use module level loggers, and always log exception tracebacks.
- Improve API documentation: separate pages per modules, each with an inheritance diagram as an overview.

### 7.2 0.6.0 (2021-04-23)

- Technix capacitor charger using either serial connection or Telnet protocol.
- **Extensions, improvements and fixes in existing devices:**
  - In `dev.tiepie.TiePieOscilloscope`:
    - \* redesigned measurement start and data collection API, incl. time out argument, with no/infinite time out option;
    - \* trigger allows now a no/infinite time out;
    - \* record length and trigger level were fixed to accept, respectively, floating point and integer numbers;
    - \* fixed resolution validation bug;
  - `dev.heinzinger.HeinzingerDI` and `dev.rs_rto1024.RTO1024` instances are now resilient to multiple `stop()` calls.
  - In `dev.crylas.CryLasLaser`: default configuration timeout and polling period were adjusted;
  - Fixed PSI9080 example script.
- **Package and source code improvements:**

- Update to backward-incompatible `pyvisa-py>=0.5.2`. Developers, do update your local development environments!
- External libraries, like LibTiePie SDK or LJM Library, are now not installed by default; they are now extra installation options.
- Added Python 3.9 support.
- Improved number formatting in logs.
- Typing improvements and fixes for `mypy>=0.800`.

## 7.3 0.5.0 (2020-11-11)

- TiePie USB oscilloscope, generator and I2C host devices, as a wrapper of the Python bindings for the LibTiePie SDK.
- a FuG Elektronik Power Supply (e.g. Capacitor Charger HCK) using the built-in ADDAT controller with the Probus V protocol over a serial connection
- All devices polling status or measurements use now a `dev.utils.Poller` utility class.
- **Extensions and improvements in existing devices:**
  - In `dev.rs_rto1024.RTO1024`: added Channel state, scale, range, position and offset accessors, and measurements activation and read methods.
  - In `dev.sst_luminox.Luminox`: added querying for all measurements in polling mode, and made output mode activation more robust.
  - In `dev.newport.NewportSMC100PP`: an error-prone `wait_until_move_finished` method was replaced by a fixed waiting time, device operations are now robust to a power supply cut, and device restart is not required to apply a start configuration.
- **Other minor improvements:**
  - Single failure-safe starting and stopping of devices sequenced via `dev.base.DeviceSequenceMixin`.
  - Moved `read_text_nonempty` up to `comm.serial.SerialCommunication`.
  - Added development Dockerfile.
  - Updated package and development dependencies: `pymodbus`, `pytest-mock`.

## 7.4 0.4.0 (2020-07-16)

- **Significantly improved new Supercube device controller:**
  - more robust error-handling,
  - status polling with generic `Poller` helper,
  - messages and status boards.
  - tested with a physical device,
- Improved OPC UA client wrapper, with better error handling, incl. re-tries on `concurrent.futures.TimeoutError`.
- SST Luminox Oxygen sensor device controller.

- **Backward-incompatible changes:**

- `CommunicationProtocol.access_lock` has changed type from `threading.Lock` to `threading.RLock`.
- `ILS2T.relative_step` and `ILS2T.absolute_position` are now called, respectively, `ILS2T.write_relative_step` and `ILS2T.write_absolute_position`.

- **Minor bugfixes and improvements:**

- fix use of max resolution in `Labjack.set_ain_resolution()`,
- resolve ILS2T devices relative and absolute position setters race condition,
- added acoustic horn function in the 2015 Supercube.

- **Toolchain changes:**

- add Python 3.8 support,
- drop pytest-runner support,
- ensure compatibility with `labjack_ljm` 2019 version library.

## 7.5 0.3.5 (2020-02-18)

- Fix issue with reading integers from LabJack LJM Library (device's product ID, serial number etc.)
- Fix development requirements specification (tox version).

## 7.6 0.3.4 (2019-12-20)

- **New devices using serial connection:**

- Heinzinger Digital Interface I/II and a Heinzinger PNC power supply
- Q-switched Pulsed Laser and a laser attenuator from CryLas
- Newport SMC100PP single axis motion controller for 2-phase stepper motors
- Pfeiffer TPG controller (TPG 25x, TPG 26x and TPG 36x) for Compact pressure Gauges

- PEP 561 compatibility and related corrections for static type checking (now in CI)

- **Refactorings:**

- Protected non-thread safe read and write in communication protocols
- Device sequence mixin: start/stop, add/rm and lookup
- `.format()` to f-strings
- more enumerations and a quite some improvements of existing code

- Improved error docstrings (`:raises:` annotations) and extended tests for errors.

## 7.7 0.3.3 (2019-05-08)

- Use PyPI labjack-ljm (no external dependencies)

## 7.8 0.3.2 (2019-05-08)

- INSTALLATION.rst with LJMPython prerequisite info

## 7.9 0.3.1 (2019-05-02)

- readthedocs.org support

## 7.10 0.3 (2019-05-02)

- Prevent an automatic close of VISA connection when not used.
- Rhode & Schwarz RTO 1024 oscilloscope using VISA interface over `TCP::INSTR`.
- Extended tests incl. messages sent to devices.
- Added Supercube device using an OPC UA client
- Added Supercube 2015 device using an OPC UA client (for interfacing with old system version)

## 7.11 0.2.1 (2019-04-01)

- Fix issue with LJMPython not being installed automatically with setuptools.

## 7.12 0.2.0 (2019-03-31)

- LabJack LJM Library communication wrapper and LabJack device.
- Modbus TCP communication protocol.
- Schneider Electric ILS2T stepper motor drive device.
- Elektro-Automatik PSI9000 current source device and VISA communication wrapper.
- Separate configuration classes for communication protocols and devices.
- Simple experiment manager class.

## 7.13 0.1.0 (2019-02-06)

- Communication protocol base and serial communication implementation.
- Device base and MBW973 implementation.



---

**CHAPTER  
EIGHT**

---

**INDICES AND TABLES**

- genindex
- modindex
- search



## PYTHON MODULE INDEX

### h

hvl\_ccb, 168  
hvl\_ccb.comm, 29  
hvl\_ccb.comm.base, 9  
hvl\_ccb.comm.labjack\_ljm, 13  
hvl\_ccb.comm.modbus\_tcp, 16  
hvl\_ccb.comm.opc, 18  
hvl\_ccb.comm.serial, 21  
hvl\_ccb.comm.telnet, 24  
hvl\_ccb.comm.visa, 26  
hvl\_ccb.configuration, 165  
hvl\_ccb.dev, 164  
hvl\_ccb.dev.base, 69  
hvl\_ccb.dev.crylas, 71  
hvl\_ccb.dev.ea\_psi9000, 81  
hvl\_ccb.dev.fug, 85  
hvl\_ccb.dev.heinzinger, 96  
hvl\_ccb.dev.labjack, 102  
hvl\_ccb.dev.mbw973, 107  
hvl\_ccb.dev.newport, 111  
hvl\_ccb.dev.pfeiffer\_tpg, 124  
hvl\_ccb.dev.rs\_rto1024, 130  
hvl\_ccb.dev.se\_ilis2t, 137  
hvl\_ccb.dev.sst\_luminox, 144  
hvl\_ccb.dev.supercube, 53  
hvl\_ccb.dev.supercube.base, 29  
hvl\_ccb.dev.supercube.constants, 35  
hvl\_ccb.dev.supercube.typ\_a, 49  
hvl\_ccb.dev.supercube.typ\_b, 52  
hvl\_ccb.dev.supercube2015, 68  
hvl\_ccb.dev.supercube2015.base, 53  
hvl\_ccb.dev.supercube2015.constants, 59  
hvl\_ccb.dev.supercube2015.typ\_a, 66  
hvl\_ccb.dev.technix, 148  
hvl\_ccb.dev.tiepie, 153  
hvl\_ccb.dev.utils, 161  
hvl\_ccb.dev.visa, 162  
hvl\_ccb.experiment\_manager, 167  
hvl\_ccb.utils, 165  
hvl\_ccb.utils.enum, 164  
hvl\_ccb.utils.typing, 165



# INDEX

## A

A (*HeinzingerPNC.UnitCurrent attribute*), 99  
A (*SupercubeOpcEndpoint attribute*), 48, 66  
ABSOLUTE (*TiePieOscilloscopeTriggerLevelMode attribute*), 160  
ABSOLUTE\_POSITION (*ILS2T.ActionsPtp attribute*), 138  
AC (*NewportConfigCommands attribute*), 111  
AC\_DoubleStage\_150kV (*PowerSetup attribute*), 47, 64  
AC\_DoubleStage\_200kV (*PowerSetup attribute*), 47, 64  
AC\_SingleStage\_100kV (*PowerSetup attribute*), 47, 64  
AC\_SingleStage\_50kV (*PowerSetup attribute*), 47, 64  
ACA (*TiePieOscilloscopeChannelCoupling attribute*), 158  
acceleration (*NewportSMC100PPConfig attribute*), 119  
ACCESS\_ENABLE (*ILS2TRegAddr attribute*), 143  
access\_lock (*CommunicationProtocol attribute*), 12  
ACTION\_JOG\_VALUE (*ILS2T attribute*), 138  
activate\_measurements () (*RTO1024 method*), 130  
activate\_output () (*Luminox method*), 144  
activated (*BreakdownDetection attribute*), 40, 60  
ACTIVE (*CryLasLaser.AnswersStatus attribute*), 74  
active (*OpcControl attribute*), 46  
actualsetvalue () (*FuGProbusVSetRegisters property*), 93  
ACV (*TiePieOscilloscopeChannelCoupling attribute*), 159  
adc\_mode () (*FuGProbusVMonitorRegisters property*), 92  
add\_device () (*DeviceSequenceMixin method*), 69  
add\_device () (*ExperimentManager method*), 167  
ADDR\_INCORRECT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 120  
address (*NewportSMC100PPConfig attribute*), 119  
address () (*VisaCommunicationConfig property*), 27  
address () (*VisaCommunicationConfig.InterfaceType method*), 27  
ADMODE (*FuGProbusIVCommands attribute*), 90  
Alarm0 (*AlarmText attribute*), 59  
Alarm1 (*Alarms attribute*), 36  
Alarm1 (*AlarmText attribute*), 35, 59  
Alarm10 (*Alarms attribute*), 36  
Alarm10 (*AlarmText attribute*), 35, 59  
Alarm100 (*Alarms attribute*), 36  
Alarm101 (*Alarms attribute*), 36  
Alarm102 (*Alarms attribute*), 36  
Alarm103 (*Alarms attribute*), 36  
Alarm104 (*Alarms attribute*), 36  
Alarm105 (*Alarms attribute*), 36  
Alarm106 (*Alarms attribute*), 36  
Alarm107 (*Alarms attribute*), 36  
Alarm108 (*Alarms attribute*), 36  
Alarm109 (*Alarms attribute*), 36  
Alarm11 (*Alarms attribute*), 37  
Alarm11 (*AlarmText attribute*), 35, 59  
Alarm110 (*Alarms attribute*), 37  
Alarm111 (*Alarms attribute*), 37  
Alarm112 (*Alarms attribute*), 37  
Alarm113 (*Alarms attribute*), 37  
Alarm114 (*Alarms attribute*), 37  
Alarm115 (*Alarms attribute*), 37  
Alarm116 (*Alarms attribute*), 37  
Alarm117 (*Alarms attribute*), 37  
Alarm118 (*Alarms attribute*), 37  
Alarm119 (*Alarms attribute*), 37  
Alarm12 (*Alarms attribute*), 37  
Alarm12 (*AlarmText attribute*), 35, 59  
Alarm120 (*Alarms attribute*), 37  
Alarm121 (*Alarms attribute*), 37  
Alarm122 (*Alarms attribute*), 37  
Alarm123 (*Alarms attribute*), 37  
Alarm124 (*Alarms attribute*), 37  
Alarm125 (*Alarms attribute*), 37  
Alarm126 (*Alarms attribute*), 37  
Alarm127 (*Alarms attribute*), 37  
Alarm128 (*Alarms attribute*), 37  
Alarm129 (*Alarms attribute*), 37

Alarm13 (*Alarms attribute*), 37  
Alarm13 (*AlarmText attribute*), 35, 59  
Alarm130 (*Alarms attribute*), 37  
Alarm131 (*Alarms attribute*), 37  
Alarm132 (*Alarms attribute*), 37  
Alarm133 (*Alarms attribute*), 37  
Alarm134 (*Alarms attribute*), 37  
Alarm135 (*Alarms attribute*), 37  
Alarm136 (*Alarms attribute*), 37  
Alarm137 (*Alarms attribute*), 37  
Alarm138 (*Alarms attribute*), 37  
Alarm139 (*Alarms attribute*), 37  
Alarm14 (*Alarms attribute*), 37  
Alarm14 (*AlarmText attribute*), 35, 59  
Alarm140 (*Alarms attribute*), 37  
Alarm141 (*Alarms attribute*), 37  
Alarm142 (*Alarms attribute*), 38  
Alarm143 (*Alarms attribute*), 38  
Alarm144 (*Alarms attribute*), 38  
Alarm145 (*Alarms attribute*), 38  
Alarm146 (*Alarms attribute*), 38  
Alarm147 (*Alarms attribute*), 38  
Alarm148 (*Alarms attribute*), 38  
Alarm149 (*Alarms attribute*), 38  
Alarm15 (*Alarms attribute*), 38  
Alarm15 (*AlarmText attribute*), 35  
Alarm150 (*Alarms attribute*), 38  
Alarm151 (*Alarms attribute*), 38  
Alarm16 (*Alarms attribute*), 38  
Alarm16 (*AlarmText attribute*), 35  
Alarm17 (*Alarms attribute*), 38  
Alarm17 (*AlarmText attribute*), 35, 59  
Alarm18 (*Alarms attribute*), 38  
Alarm18 (*AlarmText attribute*), 35  
Alarm19 (*Alarms attribute*), 38  
Alarm19 (*AlarmText attribute*), 35, 59  
Alarm2 (*Alarms attribute*), 38  
Alarm2 (*AlarmText attribute*), 35, 59  
Alarm20 (*Alarms attribute*), 38  
Alarm20 (*AlarmText attribute*), 35, 59  
Alarm21 (*Alarms attribute*), 38  
Alarm21 (*AlarmText attribute*), 35, 59  
Alarm22 (*Alarms attribute*), 38  
Alarm22 (*AlarmText attribute*), 35, 59  
Alarm23 (*Alarms attribute*), 38  
Alarm23 (*AlarmText attribute*), 35  
Alarm24 (*Alarms attribute*), 38  
Alarm24 (*AlarmText attribute*), 35  
Alarm25 (*Alarms attribute*), 38  
Alarm25 (*AlarmText attribute*), 35  
Alarm26 (*Alarms attribute*), 38  
Alarm26 (*AlarmText attribute*), 36  
Alarm27 (*Alarms attribute*), 38  
Alarm28 (*Alarms attribute*), 38  
Alarm29 (*Alarms attribute*), 38  
Alarm3 (*Alarms attribute*), 38  
Alarm3 (*AlarmText attribute*), 36, 59  
Alarm30 (*Alarms attribute*), 38  
Alarm31 (*Alarms attribute*), 38  
Alarm32 (*Alarms attribute*), 38  
Alarm33 (*Alarms attribute*), 38  
Alarm34 (*Alarms attribute*), 38  
Alarm35 (*Alarms attribute*), 38  
Alarm36 (*Alarms attribute*), 38  
Alarm37 (*Alarms attribute*), 38  
Alarm38 (*Alarms attribute*), 38  
Alarm39 (*Alarms attribute*), 39  
Alarm4 (*Alarms attribute*), 39  
Alarm4 (*AlarmText attribute*), 36, 59  
Alarm40 (*Alarms attribute*), 39  
Alarm41 (*Alarms attribute*), 39  
Alarm41 (*AlarmText attribute*), 36  
Alarm42 (*Alarms attribute*), 39  
Alarm42 (*AlarmText attribute*), 36  
Alarm43 (*Alarms attribute*), 39  
Alarm43 (*AlarmText attribute*), 36  
Alarm44 (*Alarms attribute*), 39  
Alarm44 (*AlarmText attribute*), 36  
Alarm45 (*Alarms attribute*), 39  
Alarm45 (*AlarmText attribute*), 36  
Alarm46 (*Alarms attribute*), 39  
Alarm46 (*AlarmText attribute*), 36  
Alarm47 (*Alarms attribute*), 39  
Alarm47 (*AlarmText attribute*), 36  
Alarm48 (*Alarms attribute*), 39  
Alarm48 (*AlarmText attribute*), 36  
Alarm49 (*Alarms attribute*), 39  
Alarm5 (*Alarms attribute*), 39  
Alarm5 (*AlarmText attribute*), 36, 59  
Alarm50 (*Alarms attribute*), 39  
Alarm51 (*Alarms attribute*), 39  
Alarm52 (*Alarms attribute*), 39  
Alarm53 (*Alarms attribute*), 39  
Alarm54 (*Alarms attribute*), 39  
Alarm55 (*Alarms attribute*), 39  
Alarm56 (*Alarms attribute*), 39  
Alarm57 (*Alarms attribute*), 39  
Alarm58 (*Alarms attribute*), 39  
Alarm59 (*Alarms attribute*), 39  
Alarm6 (*Alarms attribute*), 39  
Alarm6 (*AlarmText attribute*), 36, 59  
Alarm60 (*Alarms attribute*), 39  
Alarm61 (*Alarms attribute*), 39  
Alarm62 (*Alarms attribute*), 39  
Alarm63 (*Alarms attribute*), 39  
Alarm64 (*Alarms attribute*), 39  
Alarm65 (*Alarms attribute*), 39  
Alarm66 (*Alarms attribute*), 39

Alarm67 (*Alarms attribute*), 39  
 Alarm68 (*Alarms attribute*), 39  
 Alarm69 (*Alarms attribute*), 39  
 Alarm7 (*Alarms attribute*), 39  
 Alarm7 (*AlarmText attribute*), 36, 59  
 Alarm70 (*Alarms attribute*), 39  
 Alarm71 (*Alarms attribute*), 40  
 Alarm72 (*Alarms attribute*), 40  
 Alarm73 (*Alarms attribute*), 40  
 Alarm74 (*Alarms attribute*), 40  
 Alarm75 (*Alarms attribute*), 40  
 Alarm76 (*Alarms attribute*), 40  
 Alarm77 (*Alarms attribute*), 40  
 Alarm78 (*Alarms attribute*), 40  
 Alarm79 (*Alarms attribute*), 40  
 Alarm8 (*Alarms attribute*), 40  
 Alarm8 (*AlarmText attribute*), 36, 59  
 Alarm80 (*Alarms attribute*), 40  
 Alarm81 (*Alarms attribute*), 40  
 Alarm82 (*Alarms attribute*), 40  
 Alarm83 (*Alarms attribute*), 40  
 Alarm84 (*Alarms attribute*), 40  
 Alarm85 (*Alarms attribute*), 40  
 Alarm86 (*Alarms attribute*), 40  
 Alarm87 (*Alarms attribute*), 40  
 Alarm88 (*Alarms attribute*), 40  
 Alarm89 (*Alarms attribute*), 40  
 Alarm9 (*Alarms attribute*), 40  
 Alarm9 (*AlarmText attribute*), 36, 59  
 Alarm90 (*Alarms attribute*), 40  
 Alarm91 (*Alarms attribute*), 40  
 Alarm92 (*Alarms attribute*), 40  
 Alarm93 (*Alarms attribute*), 40  
 Alarm94 (*Alarms attribute*), 40  
 Alarm95 (*Alarms attribute*), 40  
 Alarm96 (*Alarms attribute*), 40  
 Alarm97 (*Alarms attribute*), 40  
 Alarm98 (*Alarms attribute*), 40  
 Alarm99 (*Alarms attribute*), 40  
*Alarms* (*class in hvl\_ccb.dev.supercube.constants*), 36  
*AlarmText* (*class in hvl\_ccb.dev.supercube.constants*), 35  
*AlarmText* (*class in hvl\_ccb.dev.supercube2015.constants*), 59  
*ALL* (*TiePieOscilloscopeAutoResolutionModes attribute*), 157  
*all\_measurements* (*LuminoxMeasurementType attribute*), 146  
*all\_measurements\_types()* (*LuminoxMeasurementType class method*), 146  
*amplitude()* (*TiePieGeneratorConfig property*), 155  
*analog\_control()* (*FuGProbusVDIRegisters property*), 92  
*ANY* (*LabJack.DeviceType attribute*), 103  
*ANY* (*LJMCommunicationConfig.ConnectionType attribute*), 14  
*ANY* (*LJMCommunicationConfig.DeviceType attribute*), 14  
*ANY* (*TiePieOscilloscopeTriggerKind attribute*), 160  
*ARBITRARY* (*TiePieGeneratorSignalType attribute*), 155  
*AsyncCommunicationProtocol* (*class in hvl\_ccb.comm.base*), 9  
*AsyncCommunicationProtocolConfig* (*class in hvl\_ccb.comm.base*), 11  
*attenuation()* (*CryLasAttenuator property*), 72  
*auto* (*EarthingStickOperatingStatus attribute*), 43  
*AUTO* (*RTO1024.TriggerModes attribute*), 130  
*auto\_laser\_on* (*CryLasLaserConfig attribute*), 78  
*auto\_resolution\_mode()* (*TiePieOscilloscopeConfig property*), 159  
*AutoNumberNameEnum* (*class in hvl\_ccb.utils.enum*), 164

## B

*B* (*SupercubeOpcEndpoint attribute*), 48, 66  
*BA* (*NewportConfigCommands attribute*), 111  
*backlash\_compensation* (*NewportSMC100PPConfig attribute*), 119  
*backup\_waveform()* (*RTO1024 method*), 131  
*bar* (*PfeifferTPG.PressureUnits attribute*), 124  
*barometric\_pressure* (*LuminoxMeasurementType attribute*), 146  
*base\_velocity* (*NewportSMC100PPConfig attribute*), 119  
*baudrate* (*CryLasAttenuatorSerialCommunicationConfig attribute*), 73  
*baudrate* (*CryLasLaserSerialCommunicationConfig attribute*), 80  
*baudrate* (*FuGSerialCommunicationConfig attribute*), 95  
*baudrate* (*HeinzingerSerialCommunicationConfig attribute*), 101  
*baudrate* (*LuminoxSerialCommunicationConfig attribute*), 147  
*baudrate* (*MBW973SerialCommunicationConfig attribute*), 110  
*baudrate* (*NewportSMC100PPSerialCommunicationConfig attribute*), 122  
*baudrate* (*PfeifferTPGSerialCommunicationConfig attribute*), 129  
*baudrate* (*SerialCommunicationConfig attribute*), 22  
*BH* (*NewportConfigCommands attribute*), 111  
*board* (*VisaCommunicationConfig attribute*), 27  
*BreakdownDetection* (*class in hvl\_ccb.dev.supercube.constants*), 40  
*BreakdownDetection* (*class in hvl\_ccb.dev.supercube2015.constants*), 60

```

bytesize (CryLasAttenuatorSerialCommunication-  

Config attribute), 73
bytesize (CryLasLaserSerialCommunicationConfig  

attribute), 80
bytesize (FuGSerialCommunicationConfig attribute),  

    95
bytesize (HeinzingerSerialCommunicationConfig at-  

tribute), 101
bytesize (LuminoxSerialCommunicationConfig at-  

tribute), 147
bytesize (MBW973SerialCommunicationConfig at-  

tribute), 110
bytesize (NewportSMC100PPSerialCommunicationConfig  

attribute), 123
bytesize (PfeifferTPGSerialCommunicationConfig at-  

tribute), 129
Bytesize (SerialCommunicationConfig attribute), 22
bytesize (SerialCommunicationConfig attribute), 22

C
c (LabJack.TemperatureUnit attribute), 104
c (LabJack.ThermocoupleType attribute), 104
calibration_factor (CryLasLaserConfig at-  

tribute), 78
calibration_mode () (FuGProbusVDIRegisters  

property), 92
cc_mode () (FuGProbusVDIRegisters property), 92
cee16 (GeneralSockets attribute), 44, 62
channels_enabled () (TiePieOscilloscope prop-  

erty), 156
check_for_error () (New-  

portSMC100PPSerialCommunication method),  

    120
check_generic_type () (in module  

hvl_ccb.utils.typing), 165
check_master_slave_config () (PSI9000  

method), 81
chunk_size (VisaCommunicationConfig attribute), 27
clean_amplitude () (TiePieGeneratorConfig  

method), 155
clean_auto_resolution_mode () (TiePieOscil-  

loscopeConfig static method), 159
clean_coupling () (TiePieOscilloscopeChannel-  

Config static method), 158
clean_enabled () (TiePieGeneratorConfig static  

method), 155
clean_enabled () (TiePieOscilloscopeChannelCon-  

fig static method), 158
clean_frequency () (TiePieGeneratorConfig  

method), 155
clean_input_range () (TiePieOscilloscopeChan-  

nelConfig method), 158
clean_offset () (TiePieGeneratorConfig method),  

    155
clean_clean_pre_sample_ratio () (TiePieOscillo-  

scopeConfig method), 159
clean_probe_offset () (TiePieOscilloscopeChan-  

nelConfig method), 158
clean_record_length () (TiePieOscilloscopeCon-  

fig method), 159
clean_resolution () (TiePieOscilloscopeConfig  

static method), 159
clean_safe_ground_enabled () (TiePieOscillo-  

scopeChannelConfig static method), 158
clean_sample_frequency () (TiePieOscillo-  

scopeConfig method), 159
clean_clean_signal_type () (TiePieGeneratorConfig  

static method), 155
clean_trigger_enabled () (TiePieOscillo-  

scopeChannelConfig static method), 158
clean_trigger_hysteresis () (TiePieOscillo-  

scopeChannelConfig method), 158
clean_trigger_kind () (TiePieOscilloscopeChan-  

nelConfig static method), 158
clean_trigger_level () (TiePieOscillo-  

scopeChannelConfig method), 158
clean_trigger_level_mode () (TiePieOscillo-  

scopeChannelConfig static method), 158
clean_trigger_timeout () (TiePieOscillo-  

scopeConfig method), 159
clean_values () (AsyncCommunicationProtocol-  

Config method), 11
clean_values () (CryLasAttenuatorConfig method),  

    72
clean_values () (CryLasLaserConfig method), 78
clean_values () (EmptyConfig method), 70, 166
clean_values () (FuGConfig method), 87
clean_values () (HeinzingerConfig method), 97
clean_values () (ILS2TConfig method), 141
clean_values () (LJMCommunicationConfig  

method), 15
clean_values () (LuminoxConfig method), 145
clean_values () (MBW973Config method), 109
clean_values () (ModbusTcpCommunicationConfig  

method), 17
clean_values () (NewportSMC100PPConfig  

method), 119
clean_values () (OpcUaCommunicationConfig  

method), 19
clean_values () (PfeifferTPGConfig method), 127
clean_values () (PSI9000Config method), 83
clean_values () (SerialCommunicationConfig  

method), 22
clean_values () (SupercubeConfiguration method),  

    33, 57
clean_values () (TechnixConfig method), 150
clean_values () (TelnetCommunicationConfig  

method), 25

```

clean\_values() (*TiePieDeviceConfig* method), 154  
 clean\_values() (*VisaCommunicationConfig* method), 28  
 close (*EarthingStickOperation* attribute), 43  
 close() (*CommunicationProtocol* method), 12  
 close() (*LJMCommunication* method), 13  
 close() (*ModbusTcpCommunication* method), 16  
 close() (*NullCommunicationProtocol* method), 12  
 close() (*OpcUaCommunication* method), 18  
 close() (*SerialCommunication* method), 21  
 close() (*TelnetCommunication* method), 24  
 close() (*VisaCommunication* method), 26  
 close\_shutter() (*CryLasLaser* method), 75  
 CLOSED (*CryLasLaser.AnswersShutter* attribute), 74  
 CLOSED (*CryLasLaser.ShutterStatus* attribute), 81  
 closed (*DoorStatus* attribute), 41, 60  
 closed (*EarthingStickStatus* attribute), 43, 61  
 CMD\_EXEC\_ERROR  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_CC  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_CONFIGURATION  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_DISABLE  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_HOMING  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_MOVING  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_NOT\_REFERENCED  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMD\_NOT\_ALLOWED\_PP  
     (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 120  
 CMD\_NOT\_ALLOWED\_READY  
     (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 CMR (*PfeifferTPG.SensorTypes* attribute), 125  
 CODE\_OR\_ADDR\_INVALID  
     (*NewportSMC100PPSerialCommunication.ControllerErrors* attribute), 120  
 collect\_measurement\_data() (*TiePieOscilloscope* method), 156  
 com() (*SingleCommDevice* property), 71  
 COM\_TIMEOUT (*NewportSMC100PPSerialCommunication.ControllerError* attribute), 120  
 command() (*FuGProbusIV* method), 90  
 command() (*LuminoxMeasurementType* property), 146  
 communication\_channel (*TechnixConfig* attribute), 150  
 CommunicationProtocol  
     (class in *hvl\_ccb.comm.base*), 12  
 CONFIG (*FuGProbusVRegisterGroups* attribute), 93  
 CONFIG (*NewportSMC100PP.StateMessages* attribute), 112  
 CONFIG (*NewportStates* attribute), 123  
 config() (*ConfigurationMixin* property), 165  
 config\_cls() (*AsyncCommunicationProtocol* static method), 9  
 config\_cls() (*ConfigurationMixin* static method), 166  
 config\_cls() (*CryLasAttenuator* static method), 72  
 config\_cls() (*CryLasAttenuatorSerialCommunication* static method), 73  
 config\_cls() (*CryLasLaser* static method), 75  
 config\_cls() (*CryLasLaserSerialCommunication* static method), 79  
 config\_cls() (*Device* static method), 69  
 config\_cls() (*FuGProbusIV* static method), 90  
 config\_cls() (*FuGSerialCommunication* static method), 94  
 config\_cls() (*HeinzingerDI* static method), 97  
 config\_cls() (*HeinzingerSerialCommunication* static method), 100  
 config\_cls() (*ILS2T* static method), 138  
 config\_cls() (*ILS2TModbusTcpCommunication* static method), 142  
 config\_cls() (*LJMCommunication* static method), 13  
 config\_cls() (*Luminox* static method), 144  
 config\_cls() (*LuminoxSerialCommunication* static method), 147  
 config\_cls() (*MBW973* static method), 107  
 config\_cls() (*MBW973SerialCommunication* static method), 109  
 config\_cls() (*ModbusTcpCommunication* static method), 16  
 config\_cls() (*NewportSMC100PP* static method), 113  
 config\_cls() (*NullCommunicationProtocol* static method), 12  
 config\_cls() (*OpcUaCommunication* static method), 18  
 config\_cls() (*PfeifferTPG* static method), 125  
 config\_cls() (*PfeifferTPGSerialCommunication*

*static method), 128*  
*config\_cls () (PSI9000 static method), 81*  
*config\_cls () (PSI9000VisaCommunication static method), 84*  
*config\_cls () (RTO1024 static method), 131*  
*config\_cls () (RTO1024VisaCommunication static method), 136*  
*config\_cls () (SerialCommunication static method), 21*  
*config\_cls () (Supercube2015Base static method), 54*  
*config\_cls () (SupercubeAOpcUaCommunication static method), 49, 67*  
*config\_cls () (SupercubeBase static method), 29*  
*config\_cls () (SupercubeBOpcUaCommunication static method), 52*  
*config\_cls () (SupercubeOpcUaCommunication static method), 34, 57*  
*config\_cls () (SyncCommunicationProtocol static method), 12*  
*config\_cls () (Technix static method), 148*  
*config\_cls () (TechnixSerialCommunication static method), 151*  
*config\_cls () (TechnixTelnetCommunication static method), 152*  
*config\_cls () (TelnetCommunication static method), 24*  
*config\_cls () (TiePieOscilloscope static method), 156*  
*config\_cls () (VisaCommunication static method), 26*  
*config\_cls () (VisaDevice static method), 162*  
*config\_gen (TiePieGeneratorMixin attribute), 155*  
*config\_gen (TiePieHS5 attribute), 156*  
*config\_i2c (TiePieI2CHostMixin attribute), 156*  
*config\_i2c (TiePieWS5 attribute), 160*  
*config\_osc (TiePieOscilloscope attribute), 157*  
*config\_osc\_channel\_dict (TiePieOscilloscope attribute), 157*  
*config\_status () (FuG property), 86*  
*configdataclass () (in module hvl\_ccb.configuration), 166*  
*configuration\_save\_json () (Configuration-Mixin method), 166*  
*ConfigurationMixin (class in hvl\_ccb.configuration), 165*  
*connection\_type (LJMCommunicationConfig attribute), 15*  
*contact\_range () (GeneralSupport class method), 44*  
*coupling () (TiePieOscilloscopeChannelConfig property), 158*  
*CR (FuGTerminators attribute), 96*  
*create\_serial\_port () (SerialCommunication Config method), 22*  
*create\_telnet () (TelnetCommunicationConfig method), 25*  
*CRLF (FuGTerminators attribute), 96*  
*CryLasAttenuator (class in hvl\_ccb.dev.crylas), 72*  
*CryLasAttenuatorConfig (class in hvl\_ccb.dev.crylas), 72*  
*CryLasAttenuatorError, 73*  
*CryLasAttenuatorSerialCommunication (class in hvl\_ccb.dev.crylas), 73*  
*CryLasAttenuatorSerialCommunicationConfig (class in hvl\_ccb.dev.crylas), 73*  
*CryLasLaser (class in hvl\_ccb.dev.crylas), 74*  
*CryLasLaser.AnswersShutter (class in hvl\_ccb.dev.crylas), 74*  
*CryLasLaser.AnswersStatus (class in hvl\_ccb.dev.crylas), 74*  
*CryLasLaser.LaserStatus (class in hvl\_ccb.dev.crylas), 75*  
*CryLasLaser.RepetitionRates (class in hvl\_ccb.dev.crylas), 75*  
*CryLasLaserConfig (class in hvl\_ccb.dev.crylas), 77*  
*CryLasLaserError, 78*  
*CryLasLaserNotReadyError, 78*  
*CryLasLaserPoller (class in hvl\_ccb.dev.crylas), 78*  
*CryLasLaserSerialCommunication (class in hvl\_ccb.dev.crylas), 79*  
*CryLasLaserSerialCommunicationConfig (class in hvl\_ccb.dev.crylas), 79*  
*CryLasLaserShutterStatus (class in hvl\_ccb.dev.crylas), 81*  
*CURRENT (FuGProbusIVCommands attribute), 90*  
*CURRENT (FuGReadbackChannels attribute), 94*  
*current () (FuG property), 86*  
*current () (Technix property), 148*  
*current\_lower\_limit (PSI9000Config attribute), 83*  
*current\_monitor () (FuG property), 86*  
*current\_primary (Power attribute), 47, 64*  
*current\_upper\_limit (PSI9000Config attribute), 83*  
*CurrentPosition (New-portSMC100PPConfig.HomeSearch attribute), 118*  
*cv\_mode () (FuGProbusVDIRegisters property), 92*

## D

*datachange\_notification () (OpcUaSubHandler method), 20*  
*datachange\_notification () (SupercubeSubscriptionHandler method), 35, 58*  
*date\_of\_manufacture (LuminoxMeasurementType attribute), 146*  
*DC (TiePieGeneratorSignalType attribute), 155*

DC\_DoubleStage\_280kV (*PowerSetup attribute*), 47, 64  
 DC\_SingleStage\_140kV (*PowerSetup attribute*), 47, 65  
 DC\_VOLTAGE\_TOO\_LOW (New-  
     *portSMC100PP.MotorErrors attribute*), 112  
 DCA (*TiePieOscilloscopeChannelCoupling attribute*), 159  
 DCV (*TiePieOscilloscopeChannelCoupling attribute*), 159  
 default\_com\_cls () (*CryLasAttenuator static method*), 72  
 default\_com\_cls () (*CryLasLaser static method*), 75  
 default\_com\_cls () (*FuGProbusIV static method*), 90  
 default\_com\_cls () (*HeinzingerDI static method*), 97  
 default\_com\_cls () (*ILS2T static method*), 139  
 default\_com\_cls () (*LabJack static method*), 104  
 default\_com\_cls () (*Luminox static method*), 144  
 default\_com\_cls () (*MBW973 static method*), 107  
 default\_com\_cls () (*NewportSMC100PP static method*), 113  
 default\_com\_cls () (*PfeifferTPG static method*), 125  
 default\_com\_cls () (*PSI9000 static method*), 82  
 default\_com\_cls () (*RTO1024 static method*), 131  
 default\_com\_cls () (*SingleCommDevice static method*), 71  
 default\_com\_cls () (*Supercube2015Base static method*), 54  
 default\_com\_cls () (*Supercube2015WithFU static method*), 66  
 default\_com\_cls () (*SupercubeB static method*), 52  
 default\_com\_cls () (*SupercubeBase static method*), 29  
 default\_com\_cls () (*SupercubeWithFU static method*), 50  
 default\_com\_cls () (*Technix method*), 149  
 default\_com\_cls () (*TiePieOscilloscope static method*), 157  
 default\_com\_cls () (*VisaDevice static method*), 162  
 DEFAULT\_IO\_SCANNING\_CONTROL\_VALUES  
     (*ILS2T attribute*), 138  
 default\_n\_attempts\_read\_text\_nonempty  
     (*AsyncCommunicationProtocolConfig attribute*), 11  
 default\_n\_attempts\_read\_text\_nonempty  
     (*FuGSerialCommunicationConfig attribute*), 95  
 default\_n\_attempts\_read\_text\_nonempty  
     (*HeinzingerSerialCommunicationConfig attribute*), 101  
 default\_number\_of\_recordings (*HeinzingerConfig attribute*), 97  
 Device (*class in hvl\_ccb.dev.base*), 69  
 device\_type (*LJMCommunicationConfig attribute*), 15  
 DeviceExistingException, 69  
 DeviceFailuresException, 69  
 devices\_failed\_start (*DeviceSequenceMixin attribute*), 70  
 devices\_failed\_start (*ExperimentManager attribute*), 167  
 devices\_failed\_stop (*DeviceSequenceMixin attribute*), 70  
 devices\_failed\_stop (*ExperimentManager attribute*), 167  
 DeviceSequenceMixin (*class in hvl\_ccb.dev.base*), 69  
 di () (*FuG property*), 86  
 digital\_control () (*FuGProbusVDIRegisters property*), 92  
 DIOChannel (*LabJack attribute*), 103  
 DISABLE (*NewportStates attribute*), 123  
 disable () (*ILS2T method*), 139  
 DISABLE\_FROM\_JOGGING  
     (*portSMC100PP.StateMessages attribute*), 112  
 DISABLE\_FROM\_MOVING  
     (*portSMC100PP.StateMessages attribute*), 112  
 DISABLE\_FROM\_READY  
     (*portSMC100PP.StateMessages attribute*), 112  
 DISABLED (*TiePieOscilloscopeAutoResolutionModes attribute*), 157  
 DisableEspStageCheck  
     (*portSMC100PPConfig.EspStageConfig attribute*), 118  
 DISPLACEMENT\_OUT\_OF\_LIMIT  
     (*portSMC100PPSerialCommunication.ControllerErrors attribute*), 120  
 display\_message\_board () (*SupercubeBase method*), 29  
 display\_status\_board () (*SupercubeBase method*), 29  
 do\_ioscanning\_write () (*ILS2T method*), 139  
 Door (*class in hvl\_ccb.dev.supercube.constants*), 41  
 DoorStatus  
     (*class in hvl\_ccb.dev.supercube.constants*), 41  
 DoorStatus  
     (*class in hvl\_ccb.dev.supercube2015.constants*), 60

## E

E (*LabJack.ThermocoupleType attribute*), 104  
 E0 (*FuGErrorcodes attribute*), 88  
 E1 (*FuGErrorcodes attribute*), 88  
 E10 (*FuGErrorcodes attribute*), 88  
 E100 (*FuGErrorcodes attribute*), 88  
 E106 (*FuGErrorcodes attribute*), 88  
 E11 (*FuGErrorcodes attribute*), 88  
 E115 (*FuGErrorcodes attribute*), 88  
 E12 (*FuGErrorcodes attribute*), 88  
 E125 (*FuGErrorcodes attribute*), 88  
 E13 (*FuGErrorcodes attribute*), 88  
 E135 (*FuGErrorcodes attribute*), 88  
 E14 (*FuGErrorcodes attribute*), 88  
 E145 (*FuGErrorcodes attribute*), 88  
 E15 (*FuGErrorcodes attribute*), 88  
 E155 (*FuGErrorcodes attribute*), 88  
 E16 (*FuGErrorcodes attribute*), 88  
 E165 (*FuGErrorcodes attribute*), 89  
 E2 (*FuGErrorcodes attribute*), 89  
 E206 (*FuGErrorcodes attribute*), 89  
 E306 (*FuGErrorcodes attribute*), 89  
 E4 (*FuGErrorcodes attribute*), 89  
 E5 (*FuGErrorcodes attribute*), 89  
 E504 (*FuGErrorcodes attribute*), 89  
 E505 (*FuGErrorcodes attribute*), 89  
 E6 (*FuGErrorcodes attribute*), 89  
 E666 (*FuGErrorcodes attribute*), 89  
 E7 (*FuGErrorcodes attribute*), 89  
 E8 (*FuGErrorcodes attribute*), 89  
 E9 (*FuGErrorcodes attribute*), 89  
 EarthingRod (class  
     *hvl\_ccb.dev.supercube.constants*), 41  
 EarthingRodStatus (class  
     *hvl\_ccb.dev.supercube.constants*), 41  
 EarthingStick (class  
     *hvl\_ccb.dev.supercube.constants*), 41  
 EarthingStick (class  
     *hvl\_ccb.dev.supercube2015.constants*), 60  
 EarthingStickMeta (class  
     *hvl\_ccb.dev.supercube.constants*), 43  
 EarthingStickOperatingStatus (class  
     *hvl\_ccb.dev.supercube.constants*), 43  
 EarthingStickOperation (class  
     *hvl\_ccb.dev.supercube.constants*), 43  
 EarthingStickStatus (class  
     *hvl\_ccb.dev.supercube.constants*), 43  
 EarthingStickStatus (class  
     *hvl\_ccb.dev.supercube2015.constants*), 61  
 EEPROM\_ACCESS\_ERROR (New-  
     *portSMC100PPSerialCommunication.ControllerErrors*  
     attribute), 120  
 EIGHT (*HeinzingerConfig.RecordingsEnum attribute*),  
     96  
 EIGHT\_BIT (*TiePieOscilloscopeResolution attribute*),  
     159  
 EIGHT\_HUNDRED\_MILLI\_VOLT (*TiePieOscilloscop-  
     eRange attribute*), 159  
 EIGHT\_VOLT (*TiePieOscilloscopeRange attribute*), 159  
 EIGHTBITS (*SerialCommunicationBytesize attribute*),  
     22  
 EIGHTY\_VOLT (*TiePieOscilloscopeRange attribute*),  
     159  
 EmptyConfig (class in *hvl\_ccb.configuration*), 166  
 EmptyConfig (class in *hvl\_ccb.dev.base*), 70  
 enable () (*ILS2T method*), 139  
 enabled () (*TiePieGeneratorConfig property*), 155  
 enabled () (*TiePieOscilloscopeChannelConfig prop-  
     erty*), 158  
 EnableEspStageCheck (New-  
     *portSMC100PPConfig.EspStageConfig attribute*), 118  
 encoding (AsyncCommunicationProtocolConfig  
     attribute), 11  
 encoding\_error\_handling (AsyncCommunicationProtocolConfig attribute), 11  
 EndOfRunSwitch (New-  
     *portSMC100PPConfig.HomeSearch attribute*),  
     118  
 EndOfRunSwitch\_and\_Index (New-  
     *portSMC100PPConfig.HomeSearch attribute*),  
     118  
 endpoint\_name (*OpcUaCommunicationConfig attribute*), 19  
 endpoint\_name (*SupercubeAOpcUaConfiguration at-  
     tribute*), 49, 68  
 endpoint\_name (*SupercubeBOpcUaConfiguration at-  
     tribute*), 52  
 error (*DoorStatus attribute*), 41, 60  
 in error (*EarthingStickStatus attribute*), 43, 61  
 ERROR (*ExperimentStatus attribute*), 168  
 in Error (*SafetyStatus attribute*), 48, 65  
 errorcode (*FuGError attribute*), 88  
 in Errors (class in *hvl\_ccb.dev.supercube.constants*), 44  
 Errors (class in *hvl\_ccb.dev.supercube2015.constants*),  
     62  
 ESP\_STAGE\_NAME\_INVALID (New-  
     *portSMC100PPSerialCommunication.ControllerErrors*  
     attribute), 120  
 in ETHERNET (*LJMCommunicationConfig.ConnectionType attribute*), 14  
 in EVEN (*SerialCommunicationParity attribute*), 23  
 event\_notification () (OpcUaSubHandler  
     method), 20  
 EXECUTE (*FuGProbusIVCommands attribute*), 90  
 execute\_absolute\_position () (*ILS2T method*),  
     139  
 execute\_on\_x () (*FuGProbusVConfigRegisters prop-*

erty), 91  
`execute_relative_step()` (*ILS2T method*), 139  
`EXECUTEONX` (*FuGProbusIVCommands attribute*), 90  
`exit_configuration()` (*NewportSMC100PP method*), 113  
`exit_configuration_wait_sec` (*NewportSMC100PPConfig attribute*), 119  
`experiment_blocked` (*EarthingRodStatus attribute*), 41  
`experiment_ready` (*EarthingRodStatus attribute*), 41  
`ExperimentError`, 167  
`ExperimentManager` (class *hvl\_ccb.experiment\_manager*), 167  
`ExperimentStatus` (class *hvl\_ccb.experiment\_manager*), 168  
`External` (*PowerSetup attribute*), 47, 65

**F**

`F` (*LabJack.TemperatureUnit attribute*), 104  
`failures` (*DeviceFailuresException attribute*), 69  
`FALLING` (*TiePieOscilloscopeTriggerKind attribute*), 160  
`FAST` (*ILS2T.Ref16Jog attribute*), 138  
`file_copy()` (*RTO1024 method*), 131  
`finish()` (*ExperimentManager method*), 167  
`FINISHED` (*ExperimentStatus attribute*), 168  
`FINISHING` (*ExperimentStatus attribute*), 168  
`FIRMWARE` (*FuGReadbackChannels attribute*), 94  
`FIVEBITS` (*SerialCommunicationBytesize attribute*), 22  
`FLT_INFO` (*ILS2TRegAddr attribute*), 143  
`FLT_MEM_DEL` (*ILS2TRegAddr attribute*), 143  
`FLT_MEM_RESET` (*ILS2TRegAddr attribute*), 143  
`FOLLOWING_ERROR` (*NewportSMC100PP.MotorErrors attribute*), 112  
`FOLLOWRAMP` (*FuGRampModes attribute*), 94  
`force_trigger()` (*TiePieOscilloscope method*), 157  
`force_value()` (*AsyncCommunicationProtocolConfig method*), 11  
`force_value()` (*CryLasAttenuatorConfig method*), 72  
`force_value()` (*CryLasAttenuatorSerialCommunicationConfig method*), 74  
`force_value()` (*CryLasLaserConfig method*), 78  
`force_value()` (*CryLasLaserSerialCommunicationConfig method*), 80  
`force_value()` (*EmptyConfig method*), 70, 166  
`force_value()` (*FuGConfig method*), 87  
`force_value()` (*FuGSerialCommunicationConfig method*), 95  
`force_value()` (*HeinzingerConfig method*), 97  
`force_value()` (*HeinzingerSerialCommunicationConfig method*), 101  
`force_value()` (*ILS2TConfig method*), 141

`force_value()` (*ILS2TModbusTcpCommunicationConfig method*), 142  
`force_value()` (*LJMCommunicationConfig method*), 15  
`force_value()` (*LuminoxConfig method*), 145  
`force_value()` (*LuminoxSerialCommunicationConfig method*), 147  
`force_value()` (*MBW973Config method*), 109  
`force_value()` (*MBW973SerialCommunicationConfig method*), 110  
`force_value()` (*ModbusTcpCommunicationConfig method*), 17  
`force_value()` (*NewportSMC100PPConfig method*), 119  
`force_value()` (*NewportSMC100PPSerialCommunicationConfig method*), 123  
`force_value()` (*OpcUaCommunicationConfig method*), 19  
`force_value()` (*PfeifferTPGConfig method*), 127  
`force_value()` (*PfeifferTPGSerialCommunicationConfig method*), 129  
`force_value()` (*PSI9000Config method*), 83  
`force_value()` (*PSI9000VisaCommunicationConfig method*), 84  
`force_value()` (*RTO1024Config method*), 136  
`force_value()` (*RTO1024VisaCommunicationConfig method*), 137  
`force_value()` (*SerialCommunicationConfig method*), 22  
`force_value()` (*SupercubeAOpcUaConfiguration method*), 49, 68  
`force_value()` (*SupercubeBOpcUaConfiguration method*), 52  
`force_value()` (*SupercubeConfiguration method*), 33, 57  
`force_value()` (*SupercubeOpcUaCommunicationConfig method*), 34, 57  
`force_value()` (*TechnixCommunicationConfig method*), 149  
`force_value()` (*TechnixConfig method*), 150  
`force_value()` (*TechnixSerialCommunicationConfig method*), 151  
`force_value()` (*TechnixTelnetCommunicationConfig method*), 152  
`force_value()` (*TelnetCommunicationConfig method*), 25  
`force_value()` (*TiePieDeviceConfig method*), 154  
`force_value()` (*VisaCommunicationConfig method*), 28  
`force_value()` (*VisaDeviceConfig method*), 163  
`FORTY_VOLT` (*TiePieOscilloscopeRange attribute*), 159  
`FOUR` (*HeinzingerConfig.RecordingsEnum attribute*), 96  
`FOUR_HUNDRED_MILLI_VOLT` (*TiePieOscilloscop-*

*eRange attribute), 159*

**FOUR\_VOLT** (*TiePieOscilloscopeRange attribute*), 159

**FOURTEEN\_BIT** (*TiePieOscilloscopeResolution attribute*), 160

**FREERUN** (*RTO1024.TriggerModes attribute*), 130

**frequency** (*Power attribute*), 47, 64

**frequency ()** (*TiePieGeneratorConfig property*), 155

**FRM** (*NewportConfigCommands attribute*), 111

**from\_json ()** (*ConfigurationMixin class method*), 166

**FRS** (*NewportConfigCommands attribute*), 111

**fso\_reset ()** (*Supercube2015WithFU method*), 67

**fso\_reset ()** (*SupercubeWithFU method*), 50

**FUG** (*class in hvl\_ccb.dev.fug*), 85

**FUGConfig** (*class in hvl\_ccb.dev.fug*), 87

**FugDigitalVal** (*class in hvl\_ccb.dev.fug*), 88

**FUGError**, 88

**FUGErrorcodes** (*class in hvl\_ccb.dev.fug*), 88

**FugMonitorModes** (*class in hvl\_ccb.dev.fug*), 89

**FugPolarities** (*class in hvl\_ccb.dev.fug*), 89

**FugProbusIV** (*class in hvl\_ccb.dev.fug*), 89

**FugProbusIVCommands** (*class in hvl\_ccb.dev.fug*), 90

**FugProbusV** (*class in hvl\_ccb.dev.fug*), 91

**FugProbusVConfigRegisters** (*class in hvl\_ccb.dev.fug*), 91

**FugProbusVDIRegisters** (*class in hvl\_ccb.dev.fug*), 92

**FugProbusVDORegisters** (*class in hvl\_ccb.dev.fug*), 92

**FugProbusVMonitorRegisters** (*class in hvl\_ccb.dev.fug*), 92

**FugProbusVRegisterGroups** (*class in hvl\_ccb.dev.fug*), 93

**FugProbusVSetRegisters** (*class in hvl\_ccb.dev.fug*), 93

**FugRampModes** (*class in hvl\_ccb.dev.fug*), 94

**FugReadbackChannels** (*class in hvl\_ccb.dev.fug*), 94

**FugSerialCommunication** (*class in hvl\_ccb.dev.fug*), 94

**FugSerialCommunicationConfig** (*class in hvl\_ccb.dev.fug*), 94

**FugTerminators** (*class in hvl\_ccb.dev.fug*), 96

**G**

**GeneralSockets** (*class in hvl\_ccb.dev.supercube.constants*), 44

**GeneralSockets** (*class in hvl\_ccb.dev.supercube2015.constants*), 62

**GeneralSupport** (*class in hvl\_ccb.dev.supercube.constants*), 44

**GeneralSupport** (*class in hvl\_ccb.dev.supercube2015.constants*), 62

**GeneralSupportMeta** (*class in hvl\_ccb.dev.supercube.constants*), 45

**GENERATOR** (*TiePieDeviceType attribute*), 154

**generator\_start ()** (*TiePieGeneratorMixin method*), 155

**generator\_stop ()** (*TiePieGeneratorMixin method*), 155

**GeneratorParameterLimits** (*class in hvl\_ccb.dev.tiepie*), 153

**get ()** (*AlarmText class method*), 36, 59

**get ()** (*MeasurementsDividerRatio class method*), 63

**get ()** (*MeasurementsScaledInput class method*), 64

**get\_acceleration ()** (*NewportSMC100PP method*), 113

**get\_acquire\_length ()** (*RTO1024 method*), 131

**get\_ain ()** (*LabJack method*), 104

**get\_by\_p\_id ()** (*LabJack.DeviceType class method*), 103

**get\_by\_p\_id ()** (*LJMCommunicationConfig.DeviceType class method*), 15

**get\_cal\_current\_source ()** (*LabJack method*), 104

**get\_cee16\_socket ()** (*Supercube2015Base method*), 54

**get\_cee16\_socket ()** (*SupercubeBase method*), 29

**get\_channel\_offset ()** (*RTO1024 method*), 131

**get\_channel\_position ()** (*RTO1024 method*), 131

**get\_channel\_range ()** (*RTO1024 method*), 131

**get\_channel\_scale ()** (*RTO1024 method*), 131

**get\_channel\_state ()** (*RTO1024 method*), 132

**get\_controller\_information ()** (*NewportSMC100PP method*), 113

**get\_current ()** (*HeinzingerDI method*), 98

**get\_dc\_volt ()** (*ILS2T method*), 139

**get\_device ()** (*DeviceSequenceMixin method*), 70

**get\_device\_by\_serial\_number ()** (*in module hvl\_ccb.dev.tiepie*), 160

**get\_devices ()** (*DeviceSequenceMixin method*), 70

**get\_digital\_input ()** (*LabJack method*), 104

**get\_door\_status ()** (*Supercube2015Base method*), 54

**get\_door\_status ()** (*SupercubeBase method*), 29

**get\_earthing\_manual ()** (*Supercube2015Base method*), 54

**get\_earthing\_rod\_status ()** (*SupercubeBase method*), 30

**get\_earthing\_status ()** (*Supercube2015Base method*), 54

**get\_earthing\_stick\_manual ()** (*SupercubeBase method*), 30

**get\_earthing\_stick\_operating\_status ()** (*SupercubeBase method*), 30

**get\_earthing\_stick\_status ()** (*SupercubeBase*

*method), 30*  
 get\_error\_code () (*ILS2T method*), 139  
 get\_error\_queue () (*VisaDevice method*), 162  
 get\_frequency () (*Supercube2015WithFU method*),  
     67  
 get\_frequency () (*SupercubeWithFU method*), 50  
 get\_fso\_active ()                   (*Supercube2015WithFU*  
     *method*), 67  
 get\_fso\_active () (*SupercubeWithFU method*), 50  
 get\_full\_scale\_mbar () (*PfeifferTPG method*),  
     125  
 get\_full\_scale\_unitless ()        (*PfeifferTPG*  
     *method*), 125  
 get\_identification () (*VisaDevice method*), 162  
 get\_interface\_version ()         (*HeinzingerDI*  
     *method*), 98  
 get\_max\_voltage ()                (*Supercube2015WithFU*  
     *method*), 67  
 get\_max\_voltage () (*SupercubeWithFU method*),  
     50  
 get\_measurement\_ratio ()         (*Supercube2015Base*  
     *method*), 54  
 get\_measurement\_ratio ()         (*SupercubeBase*  
     *method*), 30  
 get\_measurement\_voltage ()        (*Supercube2015Base*  
     *method*), 54  
 get\_measurement\_voltage ()        (*SupercubeBase*  
     *method*), 30  
 get\_motor\_configuration ()        (*New-*  
     *portSMC100PP method*), 114  
 get\_move\_duration ()             (*NewportSMC100PP*  
     *method*), 114  
 get\_negative\_software\_limit ()    (*New-*  
     *portSMC100PP method*), 114  
 get\_number\_of\_recordings ()      (*HeinzingerDI*  
     *method*), 98  
 get\_output () (*PSI9000 method*), 82  
 get\_position () (*ILS2T method*), 139  
 get\_position () (*NewportSMC100PP method*), 114  
 get\_positive\_software\_limit ()    (*New-*  
     *portSMC100PP method*), 114  
 get\_power\_setup ()                (*Supercube2015WithFU*  
     *method*), 67  
 get\_power\_setup () (*SupercubeWithFU method*),  
     50  
 get\_primary\_current ()          (*Supercube2015WithFU*  
     *method*), 67  
 get\_primary\_current ()          (*SupercubeWithFU*  
     *method*), 50  
 get\_primary\_voltage ()         (*Supercube2015WithFU*  
     *method*), 67  
 get\_primary\_voltage ()         (*SupercubeWithFU*  
     *method*), 51  
 get\_product\_id () (*LabJack method*), 105  
 get\_product\_name () (*LabJack method*), 105  
 get\_product\_type () (*LabJack method*), 105  
 get\_pulse\_energy\_and\_rate ()    (*CryLasLaser*  
     *method*), 75  
 get\_reference\_point () (*RTO1024 method*), 132  
 get\_register () (*FuGProbusV method*), 91  
 get\_repetitions () (*RTO1024 method*), 132  
 get\_sbus\_rh () (*LabJack method*), 105  
 get\_sbus\_temp () (*LabJack method*), 105  
 get\_serial\_number () (*HeinzingerDI method*), 98  
 get\_serial\_number () (*LabJack method*), 105  
 get\_state () (*NewportSMC100PP method*), 115  
 get\_status () (*ILS2T method*), 140  
 get\_status () (*Supercube2015Base method*), 55  
 get\_status () (*SupercubeBase method*), 30  
 get\_status\_byte () (*Technix method*), 149  
 get\_support\_input ()             (*Supercube2015Base*  
     *method*), 55  
 get\_support\_input ()             (*SupercubeBase method*),  
     30  
 get\_support\_output ()            (*Supercube2015Base*  
     *method*), 55  
 get\_support\_output ()            (*SupercubeBase method*),  
     31  
 get\_system\_lock () (*PSI9000 method*), 82  
 get\_t13\_socket () (*Supercube2015Base method*),  
     55  
 get\_t13\_socket () (*SupercubeBase method*), 31  
 get\_target\_voltage ()          (*Supercube2015WithFU*  
     *method*), 67  
 get\_target\_voltage ()          (*SupercubeWithFU*  
     *method*), 51  
 get\_temperature () (*ILS2T method*), 140  
 get\_timestamps () (*RTO1024 method*), 132  
 get\_ui\_lower\_limits () (*PSI9000 method*), 82  
 get\_uip\_upper\_limits () (*PSI9000 method*), 82  
 get\_voltage () (*HeinzingerDI method*), 98  
 get\_voltage\_current\_setpoint () (*PSI9000*  
     *method*), 82  
 go\_home () (*NewportSMC100PP method*), 115  
 go\_to\_configuration ()         (*NewportSMC100PP*  
     *method*), 115  
 GreenNotReady (*SafetyStatus attribute*), 48, 65  
 GreenReady (*SafetyStatus attribute*), 48, 65

## H

HARDWARE (*CryLasLaser.RepetitionRates attribute*), 75  
 has\_safe\_ground () (*TiePieOscilloscopeChannel-*  
     *Config property*), 158  
 HEAD (*CryLasLaser.AnswersStatus attribute*), 74  
 HeinzingerConfig (*class in hvl\_ccb.dev.heinzinger*),  
     96  
 HeinzingerConfig.RecordingsEnum (*class in*  
     *hvl\_ccb.dev.heinzinger*), 96

HeinzingerDI (*class in hvl\_ccb.dev.heinzinger*), 97  
 HeinzingerDI.OutputStatus (*class in hvl\_ccb.dev.heinzinger*), 97  
 HeinzingerPNC (*class in hvl\_ccb.dev.heinzinger*), 99  
 HeinzingerPNC.UnitCurrent (*class in hvl\_ccb.dev.heinzinger*), 99  
 HeinzingerPNC.UnitVoltage (*class in hvl\_ccb.dev.heinzinger*), 99  
 HeinzingerPNCDeviceNotRecognizedException (*hvl\_ccb.comm.labjack\_ljm module*), 100  
 HeinzingerPNCError, 100  
 HeinzingerPNCMaxCurrentExceededException, 100  
 HeinzingerPNCMaxVoltageExceededException, 100  
 HeinzingerSerialCommunication (*class in hvl\_ccb.dev.heinzinger*), 100  
 HeinzingerSerialCommunicationConfig (*class in hvl\_ccb.dev.heinzinger*), 100  
 HIGH (*LabJack.DIOStatus attribute*), 103  
 high\_resolution() (*FuGProbusVSetRegisters property*), 93  
 home\_search\_polling\_interval (*NewportSMC100PPConfig attribute*), 119  
 home\_search\_timeout (*NewportSMC100PPConfig attribute*), 119  
 home\_search\_type (*NewportSMC100PPConfig attribute*), 119  
 home\_search\_velocity (*NewportSMC100PPConfig attribute*), 119  
 HOME\_STARTED (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 120  
 HomeSwitch (*NewportSMC100PPConfig.HomeSearch attribute*), 118  
 HomeSwitch\_and\_Index (*NewportSMC100PPConfig.HomeSearch attribute*), 118  
 HOMING (*NewportStates attribute*), 123  
 HOMING\_FROM\_RS232 (*NewportSMC100PP.StateMessages attribute*), 112  
 HOMING\_FROM\_SMC (*NewportSMC100PP.StateMessages attribute*), 112  
 HOMING\_TIMEOUT (*NewportSMC100PP.MotorErrors attribute*), 112  
 horn (*Safety attribute*), 65  
 horn () (*Supercube2015Base method*), 55  
 host (*ModbusTcpCommunicationConfig attribute*), 17  
 host (*OpcUaCommunicationConfig attribute*), 19  
 host (*TelnetCommunicationConfig attribute*), 25  
 host (*VisaCommunicationConfig attribute*), 28  
 hPascal (*PfeifferTPG.PressureUnits attribute*), 124  
 HT (*NewportConfigCommands attribute*), 111  
 hv () (*Technix property*), 149  
 hvl\_ccb  
     module, 168  
 hvl\_ccb.comm  
     module, 29  
 hvl\_ccb.comm.base  
     module, 9  
 hvl\_ccb.comm.labjack\_ljm  
     module, 13  
 hvl\_ccb.comm.modbus\_tcp  
 hvl\_ccb.comm.opc  
 hvl\_ccb.comm.serial  
     module, 21  
 hvl\_ccb.comm.telnet  
     module, 24  
 hvl\_ccb.comm.visa  
     module, 26  
 hvl\_ccb.configuration  
     module, 165  
 hvl\_ccb.dev  
     module, 164  
 hvl\_ccb.dev.base  
     module, 69  
 hvl\_ccb.dev.crylas  
     module, 71  
 hvl\_ccb.dev.ea\_psi9000  
     module, 81  
 hvl\_ccb.dev.fug  
     module, 85  
 hvl\_ccb.dev.heinzinger  
     module, 96  
 hvl\_ccb.dev.labjack  
     module, 102  
 hvl\_ccb.dev.mbw973  
     module, 107  
 hvl\_ccb.dev.newport  
     module, 111  
 hvl\_ccb.dev.pfeiffer\_tpg  
     module, 124  
 hvl\_ccb.dev.rs\_rt01024  
     module, 130  
 hvl\_ccb.dev.se\_ilis2t  
     module, 137  
 hvl\_ccb.dev.sst\_luminox  
     module, 144  
 hvl\_ccb.dev.supercube  
     module, 53  
 hvl\_ccb.dev.supercube.base  
     module, 29  
 hvl\_ccb.dev.supercube.constants  
     module, 35

hvl\_ccb.dev.supercube.typ\_a  
     module, 49  
 hvl\_ccb.dev.supercube.typ\_b  
     module, 52  
 hvl\_ccb.dev.supercube2015  
     module, 68  
 hvl\_ccb.dev.supercube2015.base  
     module, 53  
 hvl\_ccb.dev.supercube2015.constants  
     module, 59  
 hvl\_ccb.dev.supercube2015.typ\_a  
     module, 66  
 hvl\_ccb.dev.technix  
     module, 148  
 hvl\_ccb.dev.tiepie  
     module, 153  
 hvl\_ccb.dev.utils  
     module, 161  
 hvl\_ccb.dev.visa  
     module, 162  
 hvl\_ccb.experiment\_manager  
     module, 167  
 hvl\_ccb.utils  
     module, 165  
 hvl\_ccb.utils.enum  
     module, 164  
 hvl\_ccb.utils.typing  
     module, 165  
 hysteresis\_compensation  
     (*NewportSMC100PPConfig attribute*), 119

|

I2C (*TiePieDeviceType attribute*), 154  
 I2CHostParameterLimits  
     (class in  
         *hvl\_ccb.dev.tiepie*), 153  
 ID (*FuGProbusIVCommands attribute*), 90  
 Identification\_error (*PfeifferTPG.SensorStatus attribute*), 125  
 identifier (*LJMCommunicationConfig attribute*), 15  
 identify\_device () (*FuG method*), 86  
 identify\_device () (*HeinzingerPNC method*), 99  
 identify\_sensors () (*PfeifferTPG method*), 126  
 IKR (*PfeifferTPG.SensorTypes attribute*), 125  
 IKR11 (*PfeifferTPG.SensorTypes attribute*), 125  
 IKR9 (*PfeifferTPG.SensorTypes attribute*), 125  
 ILS2T (class in *hvl\_ccb.dev.se\_ilst*), 137  
 ILS2T.ActionsPtp (class in *hvl\_ccb.dev.se\_ilst*), 138  
 ILS2T.Mode (class in *hvl\_ccb.dev.se\_ilst*), 138  
 ILS2T.Ref16Jog (class in *hvl\_ccb.dev.se\_ilst*), 138  
 ILS2T.State (class in *hvl\_ccb.dev.se\_ilst*), 138  
 ILS2TConfig (class in *hvl\_ccb.dev.se\_ilst*), 141  
 ILS2TException, 142

ILS2TModbusTcpCommunication (class in  
     *hvl\_ccb.dev.se\_ilst*), 142  
 ILS2TModbusTcpCommunicationConfig (class  
     in *hvl\_ccb.dev.se\_ilst*), 142  
 ILS2TRegAddr (class in *hvl\_ccb.dev.se\_ilst*), 143  
 ILS2TRegDatatype (class in *hvl\_ccb.dev.se\_ilst*),  
     143  
 IMMEDIATELY (*FuGRampModes attribute*), 94  
 IMR (*PfeifferTPG.SensorTypes attribute*), 125  
 in\_1\_1 (*GeneralSupport attribute*), 44, 62  
 in\_1\_2 (*GeneralSupport attribute*), 44, 62  
 in\_2\_1 (*GeneralSupport attribute*), 44, 62  
 in\_2\_2 (*GeneralSupport attribute*), 44, 62  
 in\_3\_1 (*GeneralSupport attribute*), 44, 62  
 in\_3\_2 (*GeneralSupport attribute*), 44, 62  
 in\_4\_1 (*GeneralSupport attribute*), 44, 62  
 in\_4\_2 (*GeneralSupport attribute*), 44, 62  
 in\_5\_1 (*GeneralSupport attribute*), 44, 62  
 in\_5\_2 (*GeneralSupport attribute*), 44, 62  
 in\_6\_1 (*GeneralSupport attribute*), 45, 62  
 in\_6\_2 (*GeneralSupport attribute*), 45, 62  
 INACTIVE (*CryLasLaser.AnswersStatus attribute*), 75  
 inactive (*DoorStatus attribute*), 41, 60  
 inactive (*EarthingStickStatus attribute*), 43, 61  
 inhibit () (*Technix property*), 149  
 init\_attenuation  
     (*CryLasAttenuatorConfig attribute*), 72  
 init\_monitored\_nodes ()  
     (*OpcUaCommunication method*), 18  
 init\_shutter\_status  
     (*CryLasLaserConfig attribute*), 78  
 initialize () (*NewportSMC100PP method*), 115  
 INITIALIZED (*ExperimentStatus attribute*), 168  
 INITIALIZING (*ExperimentStatus attribute*), 168  
 Initializing (*SafetyStatus attribute*), 48, 65  
 INPUT (*FuGProbusVRegisterGroups attribute*), 93  
 input () (*GeneralSupport class method*), 45, 63  
 input\_1 (*MeasurementsDividerRatio attribute*), 45, 63  
 input\_1 (*MeasurementsScaledInput attribute*), 46, 64  
 input\_2 (*MeasurementsDividerRatio attribute*), 46  
 input\_2 (*MeasurementsScaledInput attribute*), 46, 64  
 input\_3 (*MeasurementsDividerRatio attribute*), 46  
 input\_3 (*MeasurementsScaledInput attribute*), 46, 64  
 input\_4 (*MeasurementsDividerRatio attribute*), 46  
 input\_4 (*MeasurementsScaledInput attribute*), 46, 64  
 input\_range ()  
     (*TiePieOscilloscopeChannelConfig property*), 158  
 INT32 (*ILS2TRegDatatype attribute*), 143  
 interface\_type (*PSI9000VisaCommunicationConfig attribute*), 85  
 interface\_type (*RTO1024VisaCommunicationConfig attribute*), 137  
 interface\_type  
     (*VisaCommunicationConfig attribute*), 28

internal (*LabJack.CjcType attribute*), 103  
 Internal (*PowerSetup attribute*), 47, 65  
 InvalidSupercubeStatusError, 53  
 IO\_SCANNING (*ILS2TRegAddr attribute*), 143  
 IoScanningModeValueError, 143  
 is\_configdataclass (*AsyncCommunicationProtocolConfig attribute*), 11  
 is\_configdataclass (*CryLasAttenuatorConfig attribute*), 73  
 is\_configdataclass (*CryLasLaserConfig attribute*), 78  
 is\_configdataclass (*EmptyConfig attribute*), 71, 166  
 is\_configdataclass (*FuGConfig attribute*), 87  
 is\_configdataclass (*HeinzingerConfig attribute*), 97  
 is\_configdataclass (*ILS2TConfig attribute*), 141  
 is\_configdataclass (*LJMCommunicationConfig attribute*), 15  
 is\_configdataclass (*LuminoxConfig attribute*), 145  
 is\_configdataclass (*MBW973Config attribute*), 109  
 is\_configdataclass (*ModbusTcpCommunicationConfig attribute*), 17  
 is\_configdataclass (*NewportSMC100PPConfig attribute*), 119  
 is\_configdataclass (*OpcUaCommunicationConfig attribute*), 19  
 is\_configdataclass (*PfeifferTPGConfig attribute*), 127  
 is\_configdataclass (*SupercubeConfiguration attribute*), 33, 57  
 is\_configdataclass (*TechnixConfig attribute*), 150  
 is\_configdataclass (*TiePieDeviceConfig attribute*), 154  
 is\_configdataclass (*VisaCommunicationConfig attribute*), 28  
 is\_data\_ready\_polling\_interval\_sec (*TiePieDeviceConfig attribute*), 154  
 is\_done () (*MBW973 method*), 108  
 is\_error () (*ExperimentManager method*), 168  
 is\_finished () (*ExperimentManager method*), 168  
 is\_generic\_type\_hint () (in *module hvl\_ccb.utils.typing*), 165  
 is\_in\_range () (*ILS2TRegDatatype method*), 143  
 is\_inactive () (*CryLasLaser.LaserStatus property*), 75  
 is\_measurement\_data\_ready () (*TiePieOscilloscope method*), 157  
 is\_open () (*LJMCommunication property*), 13  
 is\_open () (*OpcUaCommunication property*), 18  
 is\_open () (*SerialCommunication property*), 21  
 is\_open () (*TelnetCommunication property*), 24  
 is\_polling () (*Poller method*), 161  
 is\_ready () (*CryLasLaser.LaserStatus property*), 75  
 is\_running () (*ExperimentManager method*), 168  
 is\_triggered () (*TiePieOscilloscope method*), 157  
 is\_valid\_scale\_range\_reversed\_str () (*PfeifferTPGConfig.Model method*), 127

**J**

J (*LabJack.ThermocoupleType attribute*), 104  
 jerk\_time (*NewportSMC100PPConfig attribute*), 119  
 JOG (*ILS2T.Mode attribute*), 138  
 jog\_run () (*ILS2T method*), 140  
 jog\_stop () (*ILS2T method*), 140  
 JOGGING (*NewportStates attribute*), 123  
 JOGGING\_FROM\_DISABLE (*NewportSMC100PP.StateMessages attribute*), 112  
 JOGGING\_FROM\_READY (*NewportSMC100PP.StateMessages attribute*), 112  
 JOGN\_FAST (*ILS2TRegAddr attribute*), 143  
 JOGN\_SLOW (*ILS2TRegAddr attribute*), 143  
 JR (*NewportConfigCommands attribute*), 111

**K**

K (*LabJack.TemperatureUnit attribute*), 104  
 K (*LabJack.ThermocoupleType attribute*), 104  
 keys () (*AsyncCommunicationProtocolConfig class method*), 11  
 keys () (*CryLasAttenuatorConfig class method*), 73  
 keys () (*CryLasAttenuatorSerialCommunicationConfig class method*), 74  
 keys () (*CryLasLaserConfig class method*), 78  
 keys () (*CryLasLaserSerialCommunicationConfig class method*), 80  
 keys () (*EmptyConfig class method*), 71, 166  
 keys () (*FuGConfig class method*), 87  
 keys () (*FuGSerialCommunicationConfig class method*), 95  
 keys () (*HeinzingerConfig class method*), 97  
 keys () (*HeinzingerSerialCommunicationConfig class method*), 101  
 keys () (*ILS2TConfig class method*), 141  
 keys () (*ILS2TModbusTcpCommunicationConfig class method*), 142  
 keys () (*LJMCommunicationConfig class method*), 15  
 keys () (*LuminoxConfig class method*), 145  
 keys () (*LuminoxSerialCommunicationConfig class method*), 147  
 keys () (*MBW973Config class method*), 109  
 keys () (*MBW973SerialCommunicationConfig class method*), 110  
 keys () (*ModbusTcpCommunicationConfig class method*), 17

keys () (*NewportSMC100PPConfig class method*), 119  
 keys () (*NewportSMC100PPSerialCommunicationConfig class method*), 123  
 keys () (*OpcUaCommunicationConfig class method*), 19  
 keys () (*PfeifferTPGConfig class method*), 127  
 keys () (*PfeifferTPGSerialCommunicationConfig class method*), 129  
 keys () (*PSI9000Config class method*), 84  
 keys () (*PSI9000VisaCommunicationConfig class method*), 85  
 keys () (*RTO1024Config class method*), 136  
 keys () (*RTO1024VisaCommunicationConfig class method*), 137  
 keys () (*SerialCommunicationConfig class method*), 23  
 keys () (*SupercubeAOpcUaConfiguration method*), 50, 68  
 keys () (*SupercubeBOpcUaConfiguration method*), 53  
 keys () (*SupercubeConfiguration class method*), 33, 57  
 keys () (*SupercubeOpcUaCommunicationConfig class method*), 34, 58  
 keys () (*TechnixCommunicationConfig class method*), 149  
 keys () (*TechnixConfig class method*), 150  
 keys () (*TechnixSerialCommunicationConfig class method*), 151  
 keys () (*TechnixTelnetCommunicationConfig class method*), 152  
 keys () (*TelnetCommunicationConfig class method*), 25  
 keys () (*TiePieDeviceConfig class method*), 154  
 keys () (*VisaCommunicationConfig class method*), 28  
 keys () (*VisaDeviceConfig class method*), 163  
 kV (*HeinzingerPNC.UnitVoltage attribute*), 99

**L**

LabJack (*class in hvl\_ccb.dev.labjack*), 102  
 LabJack.AInRange (*class in hvl\_ccb.dev.labjack*), 102  
 LabJack.CalMicroAmpere (*class in hvl\_ccb.dev.labjack*), 103  
 LabJack.CjcType (*class in hvl\_ccb.dev.labjack*), 103  
 LabJack.DeviceType (*class in hvl\_ccb.dev.labjack*), 103  
 LabJack.DIOStatus (*class in hvl\_ccb.dev.labjack*), 103  
 LabJack.TemperatureUnit (*hvl\_ccb.dev.labjack*), 103  
 LabJack.ThermocoupleType (*hvl\_ccb.dev.labjack*), 104  
 LabJackError, 107  
 LabJackIdentifierDIOError, 107  
 laser\_off () (*CryLasLaser method*), 76  
 laser\_on () (*CryLasLaser method*), 76

LF (*FuGTerminators attribute*), 96  
 LFCR (*FuGTerminators attribute*), 96  
 line\_1 (*MessageBoard attribute*), 46  
 line\_10 (*MessageBoard attribute*), 46  
 line\_11 (*MessageBoard attribute*), 46  
 line\_12 (*MessageBoard attribute*), 46  
 line\_13 (*MessageBoard attribute*), 46  
 line\_14 (*MessageBoard attribute*), 46  
 line\_15 (*MessageBoard attribute*), 46  
 line\_2 (*MessageBoard attribute*), 46  
 line\_3 (*MessageBoard attribute*), 46  
 line\_4 (*MessageBoard attribute*), 46  
 line\_5 (*MessageBoard attribute*), 46  
 line\_6 (*MessageBoard attribute*), 46  
 line\_7 (*MessageBoard attribute*), 46  
 line\_8 (*MessageBoard attribute*), 46  
 line\_9 (*MessageBoard attribute*), 46  
 list\_devices () (*TiePieOscilloscope static method*), 157  
 list\_directory () (*RTO1024 method*), 132  
 live (*OpcControl attribute*), 46  
 LJMCommunication (*class in hvl\_ccb.comm.labjack\_ljm*), 13  
 LJMCommunicationConfig (*class in hvl\_ccb.comm.labjack\_ljm*), 14  
 LJMCommunicationConfig.ConnectionType (*class in hvl\_ccb.comm.labjack\_ljm*), 14  
 LJMCommunicationConfig.DeviceType (*class in hvl\_ccb.comm.labjack\_ljm*), 14  
 LJMCommunicationError, 15  
 lm34 (*LabJack.CjcType attribute*), 103  
 load\_configuration () (*RTO1024 method*), 132  
 local\_display () (*RTO1024 method*), 132  
 locked (*DoorStatus attribute*), 41, 60  
 log\_set () (*in module hvl\_ccb.dev.tiepie*), 161  
 LOW (*LabJack.DIOStatus attribute*), 103  
 Luminox (*class in hvl\_ccb.dev.sst\_luminox*), 144  
 LuminoxConfig (*class in hvl\_ccb.dev.sst\_luminox*), 145  
 LuminoxMeasurementType (*class in hvl\_ccb.dev.sst\_luminox*), 146  
 LuminoxMeasurementTypeDict (*in module hvl\_ccb.dev.sst\_luminox*), 146  
 LuminoxMeasurementTypeError, 146  
 LuminoxMeasurementTypeValue (*in module hvl\_ccb.dev.sst\_luminox*), 146  
 LuminoxOutputMode (*class in hvl\_ccb.dev.sst\_luminox*), 146  
 LuminoxOutputModeError, 147  
 LuminoxSerialCommunication (*class in hvl\_ccb.dev.sst\_luminox*), 147  
 LuminoxSerialCommunicationConfig (*class in hvl\_ccb.dev.sst\_luminox*), 147

## M

mA (*HeinzingerPNC.UnitCurrent attribute*), 99  
 manual (*EarthingStickOperatingStatus attribute*), 43  
 manual () (*EarthingStick class method*), 42, 60  
 manual\_1 (*EarthingStick attribute*), 42, 60  
 manual\_2 (*EarthingStick attribute*), 42, 60  
 manual\_3 (*EarthingStick attribute*), 42, 60  
 manual\_4 (*EarthingStick attribute*), 42, 60  
 manual\_5 (*EarthingStick attribute*), 42, 60  
 manual\_6 (*EarthingStick attribute*), 42, 60  
 manuals () (*EarthingStick class method*), 42  
 MARK (*SerialCommunicationParity attribute*), 23  
 max\_current (*TechnixConfig attribute*), 150  
 max\_current () (*FuG property*), 86  
 max\_current () (*HeinzingerPNC property*), 99  
 max\_current () (*Technix property*), 149  
 max\_current.hardware () (*FuG property*), 86  
 max\_current.hardware () (*HeinzingerPNC property*), 99  
 max\_timeout\_retry\_nr (*OpcUaCommunicationConfig attribute*), 19  
 max\_voltage (*TechnixConfig attribute*), 150  
 max\_voltage () (*FuG property*), 86  
 max\_voltage () (*HeinzingerPNC property*), 99  
 max\_voltage () (*Technix property*), 149  
 max\_voltage.hardware () (*FuG property*), 86  
 max\_voltage.hardware () (*HeinzingerPNC property*), 100  
 mbar (*PfeifferTPG.PressureUnits attribute*), 124  
 MBW973 (*class in hvl\_ccb.dev.mbw973*), 107  
 MBW973Config (*class in hvl\_ccb.dev.mbw973*), 108  
 MBW973ControlRunningException, 109  
 MBW973Error, 109  
 MBW973PumpRunningException, 109  
 MBW973SerialCommunication (*class in hvl\_ccb.dev.mbw973*), 109  
 MBW973SerialCommunicationConfig (*class in hvl\_ccb.dev.mbw973*), 109  
 measure () (*PfeifferTPG method*), 126  
 measure\_all () (*PfeifferTPG method*), 126  
 measure\_current () (*HeinzingerDI method*), 98  
 measure\_voltage () (*HeinzingerDI method*), 98  
 measure\_voltage\_current () (*PSI9000 method*), 82  
 MeasurementsDividerRatio (*class in hvl\_ccb.dev.supercube.constants*), 45  
 MeasurementsDividerRatio (*class in hvl\_ccb.dev.supercube2015.constants*), 63  
 MeasurementsScaledInput (*class in hvl\_ccb.dev.supercube.constants*), 46  
 MeasurementsScaledInput (*class in hvl\_ccb.dev.supercube2015.constants*), 63  
 message (*Errors attribute*), 44  
 MessageBoard (*class in hvl\_ccb.dev.supercube.constants*), 46  
 micro\_step\_per\_full\_step\_factor (*NewportSMC100PPConfig attribute*), 119  
 Micron (*PfeifferTPG.PressureUnits attribute*), 124  
 ModbusTcpCommunication (*class in hvl\_ccb.comm.modbus\_tcp*), 16  
 ModbusTcpCommunicationConfig (*class in hvl\_ccb.comm.modbus\_tcp*), 17  
 ModbusTcpConnectionFailedException, 17  
 model (*PfeifferTPGConfig attribute*), 127  
 module  
     hvl\_ccb, 168  
     hvl\_ccb.comm, 29  
     hvl\_ccb.comm.base, 9  
     hvl\_ccb.comm.labjack\_ljm, 13  
     hvl\_ccb.comm.modbus\_tcp, 16  
     hvl\_ccb.comm.opc, 18  
     hvl\_ccb.comm.serial, 21  
     hvl\_ccb.comm.telnet, 24  
     hvl\_ccb.comm.visa, 26  
     hvl\_ccb.configuration, 165  
     hvl\_ccb.dev, 164  
     hvl\_ccb.dev.base, 69  
     hvl\_ccb.dev.crylas, 71  
     hvl\_ccb.dev.ea\_psi9000, 81  
     hvl\_ccb.dev.fug, 85  
     hvl\_ccb.dev.heinzinger, 96  
     hvl\_ccb.dev.labjack, 102  
     hvl\_ccb.dev.mbw973, 107  
     hvl\_ccb.dev.newport, 111  
     hvl\_ccb.dev.pfeiffer\_tpg, 124  
     hvl\_ccb.dev.rs\_rto1024, 130  
     hvl\_ccb.dev.se\_ilis2t, 137  
     hvl\_ccb.dev.sst\_luminox, 144  
     hvl\_ccb.dev.supercube, 53  
     hvl\_ccb.dev.supercube.base, 29  
     hvl\_ccb.dev.supercube.constants, 35  
     hvl\_ccb.dev.supercube.typ\_a, 49  
     hvl\_ccb.dev.supercube.typ\_b, 52  
     hvl\_ccb.dev.supercube2015, 68  
     hvl\_ccb.dev.supercube2015.base, 53  
     hvl\_ccb.dev.supercube2015.constants, 59  
     hvl\_ccb.dev.supercube2015.typ\_a, 66  
     hvl\_ccb.dev.technix, 148  
     hvl\_ccb.dev.tiepie, 153  
     hvl\_ccb.dev.utils, 161  
     hvl\_ccb.dev.visa, 162  
     hvl\_ccb.experiment\_manager, 167  
     hvl\_ccb.utils, 165  
     hvl\_ccb.utils.enum, 164  
     hvl\_ccb.utils.typing, 165  
 MONITOR\_I (*FuGProbusVRegisterGroups attribute*), 93

MONITOR\_V (*FuGProbusVRegisterGroups attribute*), 93  
most\_recent\_error() (*FuGProbusVConfigRegisters property*), 91  
motion\_distance\_per\_full\_step (*NewportSMC100PPConfig attribute*), 119  
motor\_config() (*NewportSMC100PPConfig property*), 119  
move\_to\_absolute\_position () (*NewportSMC100PP method*), 115  
move\_to\_relative\_position () (*NewportSMC100PP method*), 116  
move\_wait\_sec (*NewportSMC100PPConfig attribute*), 119  
MOVING (*NewportSMC100PP.StateMessages attribute*), 112  
MOVING (*NewportStates attribute*), 123  
MS\_NOMINAL\_CURRENT (*PSI9000 attribute*), 81  
MS\_NOMINAL\_VOLTAGE (*PSI9000 attribute*), 81  
msb\_first () (*TechnixStatusByte method*), 152  
MULTI\_COMMANDS\_MAX (*VisaCommunication attribute*), 26  
MULTI\_COMMANDS\_SEPARATOR (*VisaCommunication attribute*), 26

**N**

n\_channels () (*TiePieOscilloscope property*), 157  
n\_max\_try\_get\_device (*TiePieDeviceConfig attribute*), 154  
NameEnum (*class in hvl\_ccb.utils.enum*), 164  
NAMES (*SerialCommunicationParity attribute*), 23  
names () (*RTO1024.TriggerModes class method*), 130  
namespace\_index (*SupercubeConfiguration attribute*), 33, 57  
NATIVEONLY (*TiePieOscilloscopeAutoResolutionModes attribute*), 158  
NED\_END\_OF\_TURN (*NewportSMC100PP.MotorErrors attribute*), 112  
NEG (*ILS2T.Ref16Jog attribute*), 138  
NEG\_FAST (*ILS2T.Ref16Jog attribute*), 138  
NEGATIVE (*FuGPolarities attribute*), 89  
negative\_software\_limit (*NewportSMC100PPConfig attribute*), 119  
NewportConfigCommands (*class in hvl\_ccb.dev.newport*), 111  
NewportControllerError, 111  
NewportMotorError, 111  
NewportMotorPowerSupplyWasCutError, 111  
NewportSerialCommunicationError, 123  
NewportSMC100PP (*class in hvl\_ccb.dev.newport*), 112  
NewportSMC100PP.MotorErrors (*class in hvl\_ccb.dev.newport*), 112  
NewportSMC100PP.StateMessages (*class in hvl\_ccb.dev.newport*), 112  
NewportSMC100PPConfig (*class in hvl\_ccb.dev.newport*), 118  
NewportSMC100PPConfig.EspStageConfig (*class in hvl\_ccb.dev.newport*), 118  
NewportSMC100PPConfig.HomeSearch (*class in hvl\_ccb.dev.newport*), 118  
NewportSMC100PPSerialCommunication (*class in hvl\_ccb.dev.newport*), 120  
NewportSMC100PPSerialCommunication.ControllerErrors (*class in hvl\_ccb.dev.newport*), 120  
NewportSMC100PPSerialCommunicationConfig (*class in hvl\_ccb.dev.newport*), 122  
NewportStates (*class in hvl\_ccb.dev.newport*), 123  
NewportUncertainPositionError, 124  
NO (*FuDigitalVal attribute*), 88  
NO\_ERROR (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 120  
NO\_REF (*NewportStates attribute*), 123  
NO\_REF\_ESP\_STAGE\_ERROR (*NewportSMC100PP.StateMessages attribute*), 112  
NO\_REF\_FROM\_CONFIG (*NewportSMC100PP.StateMessages attribute*), 112  
NO\_REF\_FROM\_DISABLED (*NewportSMC100PP.StateMessages attribute*), 112  
NO\_REF\_FROM\_HOMING (*NewportSMC100PP.StateMessages attribute*), 112  
NO\_REF\_FROM\_JOGGING (*NewportSMC100PP.StateMessages attribute*), 112  
NO\_REF\_FROM\_MOVING (*NewportSMC100PP.StateMessages attribute*), 112  
NO\_REF\_FROM\_READY (*NewportSMC100PP.StateMessages attribute*), 113  
NO\_REF\_FROM\_RESET (*NewportSMC100PP.StateMessages attribute*), 113  
No\_sensor (*PfeifferTPG.SensorStatus attribute*), 125  
NOISE (*TiePieGeneratorSignalType attribute*), 155  
NONE (*ILS2T.Ref16Jog attribute*), 138  
NONE (*LabJack.ThermocoupleType attribute*), 104  
None (*PfeifferTPG.SensorTypes attribute*), 125  
NONE (*SerialCommunicationParity attribute*), 23  
NoPower (*PowerSetup attribute*), 48  
NORMAL (*RTO1024.TriggerModes attribute*), 130  
noSen (*PfeifferTPG.SensorTypes attribute*), 125  
noSENSOR (*PfeifferTPG.SensorTypes attribute*), 125  
not\_defined (*AlarmText attribute*), 36, 59  
nr\_trials\_activate (*LuminoxConfig attribute*),

<p>145  <b>NullCommunicationProtocol</b> (class in <i>hvl_ccb.comm.base</i>), 12  <b>Number</b> (<i>in module hvl_ccb.utils.typing</i>), 165  <b>number()</b> (<i>EarthingStick property</i>), 42  <b>number_of_decimals</b> (<i>HeinzingerConfig attribute</i>), 97  <b>number_of_sensors()</b> (<i>PfeifferTPG property</i>), 126</p> <p><b>O</b></p> <p><b>ODD</b> (<i>SerialCommunicationParity attribute</i>), 23  <b>OFF</b> (<i>FuGDigitalVal attribute</i>), 88  <b>OFF</b> (<i>HeinzingerDI.OutputStatus attribute</i>), 97  <b>offset()</b> (<i>TiePieGeneratorConfig property</i>), 155  <b>OH</b> (<i>NewportConfigCommands attribute</i>), 111  <b>Ok</b> (<i>PfeifferTPG.SensorStatus attribute</i>), 125  <b>ON</b> (<i>FuGDigitalVal attribute</i>), 88  <b>ON</b> (<i>HeinzingerDI.OutputStatus attribute</i>), 97  <b>ON</b> (<i>ILS2T.State attribute</i>), 138  <b>on()</b> (<i>FuG property</i>), 86  <b>on()</b> (<i>FuGProbusVDIRegisters property</i>), 92  <b>ONE</b> (<i>HeinzingerConfig.RecordingsEnum attribute</i>), 96  <b>ONE</b> (<i>LabJack.AInRange attribute</i>), 102  <b>ONE</b> (<i>SerialCommunicationStopbits attribute</i>), 23  <b>ONE_HUNDREDTH</b> (<i>LabJack.AInRange attribute</i>), 102  <b>ONE_POINT_FIVE</b> (<i>SerialCommunicationStopbits attribute</i>), 24  <b>ONE_TENTH</b> (<i>LabJack.AInRange attribute</i>), 103  <b>ONLYUPWARDSOFFTOZERO</b> (<i>FuGRampModes attribute</i>), 94  <b>OpcControl</b> (class in <i>hvl_ccb.dev.supercube.constants</i>), 46  <b>OpcUaCommunication</b> (class in <i>hvl_ccb.comm.opc</i>), 18  <b>OpcUaCommunicationConfig</b> (class in <i>hvl_ccb.comm.opc</i>), 19  <b>OpcUaCommunicationIOError</b>, 20  <b>OpcUaCommunicationTimeoutError</b>, 20  <b>OpcUaSubHandler</b> (class in <i>hvl_ccb.comm.opc</i>), 20  <b>open</b> (<i>DoorStatus attribute</i>), 41, 60  <b>open</b> (<i>EarthingStickOperation attribute</i>), 43  <b>open</b> (<i>EarthingStickStatus attribute</i>), 43, 62  <b>open()</b> (<i>CommunicationProtocol method</i>), 12  <b>open()</b> (<i>LJMCommunication method</i>), 13  <b>open()</b> (<i>ModbusTcpCommunication method</i>), 16  <b>open()</b> (<i>NullCommunicationProtocol method</i>), 12  <b>open()</b> (<i>OpcUaCommunication method</i>), 18  <b>open()</b> (<i>SerialCommunication method</i>), 21  <b>open()</b> (<i>TelnetCommunication method</i>), 24  <b>open()</b> (<i>VisaCommunication method</i>), 26  <b>open_shutter()</b> (<i>CryLasLaser method</i>), 76  <b>open_timeout</b> (<i>VisaCommunicationConfig attribute</i>), 28  <b>OPENED</b> (<i>CryLasLaser.AnswersShutter attribute</i>), 74</p>	<p><b>OPENED</b> (<i>CryLasLaserShutterStatus attribute</i>), 81  <b>operate()</b> (<i>Supercube2015Base method</i>), 55  <b>operate()</b> (<i>SupercubeBase method</i>), 31  <b>operate_earthing_stick()</b> (<i>SupercubeBase method</i>), 31  <b>operating_status()</b> (<i>EarthingStick class method</i>), 42  <b>operating_status_1</b> (<i>EarthingStick attribute</i>), 42  <b>operating_status_2</b> (<i>EarthingStick attribute</i>), 42  <b>operating_status_3</b> (<i>EarthingStick attribute</i>), 42  <b>operating_status_4</b> (<i>EarthingStick attribute</i>), 42  <b>operating_status_5</b> (<i>EarthingStick attribute</i>), 42  <b>operating_status_6</b> (<i>EarthingStick attribute</i>), 42  <b>operating_statuses()</b> (<i>EarthingStick class method</i>), 42  <b>optional_defaults()</b> (<i>AsyncCommunicationProtocolConfig class method</i>), 11  <b>optional_defaults()</b> (<i>CryLasAttenuatorConfig class method</i>), 73  <b>optional_defaults()</b> (<i>CryLasAttenuatorSerialCommunicationConfig class method</i>), 74  <b>optional_defaults()</b> (<i>CryLasLaserConfig class method</i>), 78  <b>optional_defaults()</b> (<i>CryLasLaserSerialCommunicationConfig class method</i>), 80  <b>optional_defaults()</b> (<i>EmptyConfig class method</i>), 71, 166  <b>optional_defaults()</b> (<i>FuGConfig class method</i>), 87  <b>optional_defaults()</b> (<i>FuGSerialCommunicationConfig class method</i>), 95  <b>optional_defaults()</b> (<i>HeinzingerConfig class method</i>), 97  <b>optional_defaults()</b> (<i>HeinzingerSerialCommunicationConfig class method</i>), 101  <b>optional_defaults()</b> (<i>ILS2TConfig class method</i>), 141  <b>optional_defaults()</b> (<i>ILS2TModbusTcpCommunicationConfig class method</i>), 142  <b>optional_defaults()</b> (<i>LJMCommunicationConfig class method</i>), 15  <b>optional_defaults()</b> (<i>LuminoxConfig class method</i>), 145  <b>optional_defaults()</b> (<i>LuminoxSerialCommunicationConfig class method</i>), 147  <b>optional_defaults()</b> (<i>MBW973Config class method</i>), 109  <b>optional_defaults()</b> (<i>MBW973SerialCommunicationConfig class method</i>), 110  <b>optional_defaults()</b> (<i>ModbusTcpCommunicationConfig class method</i>), 17  <b>optional_defaults()</b> (<i>NewportSMC100PPConfig</i>)</p>
--	---

```

    class method), 119
optional_defaults() (New-
    portSMC100PPSerialCommunicationConfig
    class method), 123
optional_defaults() (OpcUaCommunication-
    Config class method), 19
optional_defaults() (PfeifferTPGConfig class
    method), 127
optional_defaults() (PfeifferTPGSerialCommu-
    nicationConfig class method), 129
optional_defaults() (PSI9000Config class
    method), 84
optional_defaults() (PSI9000VisaCommunicationConfig
    method), 85
optional_defaults() (RTO1024Config class
    method), 136
optional_defaults() (RTO1024VisaCommunicationConfig class
    method), 137
optional_defaults() (SerialCommunicationCon-
    fig class method), 23
optional_defaults() (SupercubeAOpcUaConfig-
    uration class method), 50, 68
optional_defaults() (SupercubeBOpcUaConfig-
    uration class method), 53
optional_defaults() (SupercubeConfiguration
    class method), 33, 57
optional_defaults() (SupercubeOpcUaCommu-
    nicationConfig class method), 34, 58
optional_defaults() (TechnixCommunication-
    Config class method), 149
optional_defaults() (TechnixConfig class
    method), 150
optional_defaults() (TechnixSerialCommunica-
    tionConfig class method), 151
optional_defaults() (TechnixTelnetCommunica-
    tionConfig class method), 152
optional_defaults() (TelnetCommunicationCon-
    fig class method), 25
optional_defaults() (TiePieDeviceConfig class
    method), 154
optional_defaults() (VisaCommunicationConfig
    class method), 28
optional_defaults() (VisaDeviceConfig class
    method), 163
OSCILLOSCOPE (TiePieDeviceType attribute), 154
OscilloscopeChannelParameterLimits (class
    in hvl_ccb.dev.tiepie), 153
OscilloscopeParameterLimits (class in
    hvl_ccb.dev.tiepie), 153
OT (NewportConfigCommands attribute), 111
out () (FuGProbusVDORegisters property), 92
out_1_1 (GeneralSupport attribute), 45, 63
out_1_2 (GeneralSupport attribute), 45, 63
out_2_1 (GeneralSupport attribute), 45, 63
out_2_2 (GeneralSupport attribute), 45, 63
out_3_1 (GeneralSupport attribute), 45, 63
out_3_2 (GeneralSupport attribute), 45, 63
out_4_1 (GeneralSupport attribute), 45, 63
out_4_2 (GeneralSupport attribute), 45, 63
out_5_1 (GeneralSupport attribute), 45, 63
out_5_2 (GeneralSupport attribute), 45, 63
out_6_1 (GeneralSupport attribute), 45, 63
out_6_2 (GeneralSupport attribute), 45, 63
OUTPUT (FuGProbusIVCommands attribute), 90
output () (GeneralSupport class method), 45, 63
output_off () (FuGProbusIV method), 90
output_off () (HeinzingerDI method), 98
output_on () (HeinzingerDI method), 98
OUTPUT_POWER_EXCEEDED (New-
    portSMC100PP.MotorErrors attribute), 112
output_status () (HeinzingerDI property), 98
OUTPUTONCMD (FuGProbusVRegisterGroups attribute), 93
OUTPUTX0 (FuGProbusVRegisterGroups attribute), 93
OUTPUTX1 (FuGProbusVRegisterGroups attribute), 93
OUTPUTX2 (FuGProbusVRegisterGroups attribute), 93
OUTPUTXCMD (FuGProbusVRegisterGroups attribute), 93
outX0 () (FuG property), 86
outX1 () (FuG property), 86
outX2 () (FuG property), 87
outXCMD () (FuG property), 87
Overrange (PfeifferTPG.SensorStatus attribute), 125

P
PARAM_MISSING_OR_INVALID (New-
    portSMC100PPSerialCommunication.ControllerErrors
    attribute), 120
parity (CryLasAttenuatorSerialCommunicationConfig
    attribute), 74
parity (CryLasLaserSerialCommunicationConfig at-
    tribute), 80
parity (FuGSerialCommunicationConfig attribute), 95
parity (HeinzingerSerialCommunicationConfig at-
    tribute), 101
parity (LuminoxSerialCommunicationConfig at-
    tribute), 147
parity (MBW973SerialCommunicationConfig at-
    tribute), 110
parity (NewportSMC100PPSerialCommunicationConfig
    attribute), 123
parity (PfeifferTPGSerialCommunicationConfig at-
    tribute), 129
Parity (SerialCommunicationConfig attribute), 22
parity (SerialCommunicationConfig attribute), 23

```

```

parse_read_measurement_value()      (Lumi-  

    nnoxMeasurementType method), 146  

partial_pressure_o2 (LuminoxMeasurementType  

    attribute), 146  

Pascal (PfeifferTPG.PressureUnits attribute), 124  

PBR (PfeifferTPG.SensorTypes attribute), 125  

PEAK_CURRENT_LIMIT                (New-  

    portSMC100PP.MotorErrors        attribute),  

    112  

peak_output_current_limit          (New-  

    portSMC100PPConfig attribute), 119  

percent_o2 (LuminoxMeasurementType attribute),  

    146  

PfeifferTPG (class in hvl_ccb.dev.pfeiffer_tpg), 124  

PfeifferTPG.PressureUnits (class in  

    hvl_ccb.dev.pfeiffer_tpg), 124  

PfeifferTPG.SensorStatus (class in  

    hvl_ccb.dev.pfeiffer_tpg), 124  

PfeifferTPG.SensorTypes (class in  

    hvl_ccb.dev.pfeiffer_tpg), 125  

PfeifferTPGConfig (class in  

    hvl_ccb.dev.pfeiffer_tpg), 127  

PfeifferTPGConfig.Model (class in  

    hvl_ccb.dev.pfeiffer_tpg), 127  

PfeifferTPGError, 128  

PfeifferTPGSerialCommunication (class in  

    hvl_ccb.dev.pfeiffer_tpg), 128  

PfeifferTPGSerialCommunicationConfig  

    (class in hvl_ccb.dev.pfeiffer_tpg), 128  

PKR (PfeifferTPG.SensorTypes attribute), 125  

POLARITY (FuGProbusIVCommands attribute), 90  

Poller (class in hvl_ccb.dev.utils), 161  

polling (LuminoxOutputMode attribute), 146  

polling_delay_sec (SupercubeConfiguration attribute), 33  

polling_interval (MBW973Config attribute), 109  

polling_interval_sec (SupercubeConfiguration attribute), 33  

polling_interval_sec (TechnixConfig attribute),  

    150  

polling_period (CryLasLaserConfig attribute), 78  

polling_timeout (CryLasLaserConfig attribute), 78  

port (ModbusTcpCommunicationConfig attribute), 17  

port (OpcUaCommunicationConfig attribute), 20  

port (SerialCommunicationConfig attribute), 23  

port (SupercubeOpcUaCommunicationConfig attribute), 58  

port (TechnixTelnetCommunicationConfig attribute),  

    152  

port (TelnetCommunicationConfig attribute), 25  

port (VisaCommunicationConfig attribute), 28  

port_range () (GeneralSupport class method), 45  

POS (ILS2T.ReflJog attribute), 138  

POS_END_OF_TURN (NewportSMC100PP.MotorErrors  

    attribute), 112  

POS_FAST (ILS2T.ReflJog attribute), 138  

POSITION (ILS2TRegAddr attribute), 143  

POSITION_OUT_OF_LIMIT              (New-  

    portSMC100PPSerialCommunication.ControllerErrors  

    attribute), 120  

POSITIVE (FuGPolarities attribute), 89  

positive_software_limit           (New-  

    portSMC100PPConfig attribute), 119  

post_force_value () (NewportSMC100PPConfig method), 119  

post_stop_pause_sec (TechnixConfig attribute),  

    150  

Power (class in hvl_ccb.dev.supercube.constants), 47  

Power (class in hvl_ccb.dev.supercube2015.constants),  

    64  

power_limit (PSI9000Config attribute), 84  

PowerSetup (class in  

    hvl_ccb.dev.supercube.constants), 47  

PowerSetup (class in  

    hvl_ccb.dev.supercube2015.constants), 64  

pre_sample_ratio () (TiePieOscilloscopeConfig property), 159  

prepare_ultra_segmentation () (RTO1024 method), 133  

probe_offset () (TiePieOscilloscopeChannelConfig property), 158  

PSI9000 (class in hvl_ccb.dev.ea_psi9000), 81  

PSI9000Config (class in hvl_ccb.dev.ea_psi9000), 83  

PSI9000Error, 84  

PSI9000VisaCommunication (class in  

    hvl_ccb.dev.ea_psi9000), 84  

PSI9000VisaCommunicationConfig (class in  

    hvl_ccb.dev.ea_psi9000), 84  

PT100 (LabJack.ThermocoupleType attribute), 104  

PT1000 (LabJack.ThermocoupleType attribute), 104  

PT500 (LabJack.ThermocoupleType attribute), 104  

PTP (ILS2T.Mode attribute), 138  

PublicPropertiesReprMixin (class in  

    hvl_ccb.dev.tiepie), 153  

PULSE (TiePieGeneratorSignalType attribute), 155

```

## Q

```

QIL (NewportConfigCommands attribute), 111  

QUERY (FuGProbusIVCommands attribute), 90  

query () (CryLasLaserSerialCommunication method),  

    79  

query () (FuGSerialCommunication method), 94  

query () (NewportSMC100PPSerialCommunication method), 121  

query () (PfeifferTPGSerialCommunication method),  

    128  

query () (SyncCommunicationProtocol method), 12  

query () (TechnixCommunication method), 149

```

query () (*VisaCommunication method*), 27  
 query\_all () (*CryLasLaserSerialCommunication method*), 79  
 query\_multiple () (*NewportSMC100PPSerialCommunication method*), 121  
 query\_polling () (*Luminox method*), 144  
 QUICKSTOP (*ILS2T.State attribute*), 138  
 QuickStop (*SafetyStatus attribute*), 48, 65  
 quickstop () (*ILS2T method*), 140  
 quit (*Errors attribute*), 44, 62  
 quit\_error () (*Supercube2015Base method*), 55  
 quit\_error () (*SupercubeBase method*), 31

## R

R (*LabJack.ThermocoupleType attribute*), 104  
 raise\_ () (*FuGErrorcodes method*), 89  
 RAMP\_ACC (*ILS2TRegAddr attribute*), 143  
 RAMP\_DECEL (*ILS2TRegAddr attribute*), 143  
 RAMP\_N\_MAX (*ILS2TRegAddr attribute*), 143  
 RAMP\_TYPE (*ILS2TRegAddr attribute*), 143  
 rampmode () (*FuGProbusVSetRegisters property*), 93  
 ramprate () (*FuGProbusVSetRegisters property*), 93  
 rampstate () (*FuGProbusVSetRegisters property*), 93  
 RAMPUPWARDS (*FuGRampModes attribute*), 94  
 range () (*EarthingStick class method*), 42  
 RATEDCURRENT (*FuGReadbackChannels attribute*), 94  
 RATEDVOLTAGE (*FuGReadbackChannels attribute*), 94  
 read () (*AsyncCommunicationProtocol method*), 9  
 read () (*CryLasLaserSerialCommunication method*), 79  
 read () (*MBW973 method*), 108  
 read () (*OpcUaCommunication method*), 18  
 read () (*Supercube2015Base method*), 55  
 read () (*SupercubeBase method*), 31  
 read\_all () (*AsyncCommunicationProtocol method*), 10  
 read\_bytes () (*AsyncCommunicationProtocol method*), 10  
 read\_bytes () (*SerialCommunication method*), 21  
 read\_bytes () (*TelnetCommunication method*), 24  
 read\_float () (*MBW973 method*), 108  
 read\_holding\_registers () (*ModbusTcpCommunication method*), 16  
 read\_input\_registers () (*ModbusTcpCommunication method*), 16  
 read\_int () (*MBW973 method*), 108  
 read\_measurement () (*RTO1024 method*), 133  
 read\_measurements () (*MBW973 method*), 108  
 read\_name () (*LJMCommunication method*), 13  
 read\_nonempty () (*AsyncCommunicationProtocol method*), 10  
 read\_resistance () (*LabJack method*), 105

read\_single\_bytes () (*SerialCommunication method*), 21  
 read\_streaming () (*Luminox method*), 145  
 read\_termination (*VisaCommunicationConfig attribute*), 28  
 read\_text () (*AsyncCommunicationProtocol method*), 10  
 read\_text () (*NewportSMC100PPSerialCommunication method*), 121  
 read\_text\_nonempty () (*AsyncCommunicationProtocol method*), 10  
 READ\_TEXT\_SKIP\_PREFIXES (*CryLasLaserSerialCommunication attribute*), 79  
 read\_thermocouple () (*LabJack method*), 105  
 readback\_data () (*FuGProbusVConfigRegisters property*), 91  
 READBACKCHANNEL (*FuGProbusIVCommands attribute*), 90  
 READY (*CryLasLaser.AnswersStatus attribute*), 75  
 READY (*ILS2T.State attribute*), 138  
 READY (*NewportStates attribute*), 123  
 ready () (*Supercube2015Base method*), 55  
 ready () (*SupercubeBase method*), 31  
 READY\_ACTIVE (*CryLasLaser.LaserStatus attribute*), 75  
 READY\_FROM\_DISABLE (*NewportSMC100PP.StateMessages attribute*), 113  
 READY\_FROM\_HOMING (*NewportSMC100PP.StateMessages attribute*), 113  
 READY\_FROM\_JOGGING (*NewportSMC100PP.StateMessages attribute*), 113  
 READY\_FROM\_MOVING (*NewportSMC100PP.StateMessages attribute*), 113  
 READY\_INACTIVE (*CryLasLaser.LaserStatus attribute*), 75  
 record\_length () (*TiePieOscilloscopeConfig property*), 159  
 RedOperate (*SafetyStatus attribute*), 48, 66  
 RedReady (*SafetyStatus attribute*), 48, 66  
 reg\_3 () (*FuGProbusVDIRegisters property*), 92  
 RegAddr (*ILS2T attribute*), 138  
 RegDatatype (*ILS2T attribute*), 138  
 register\_pulse\_time (*TechnixConfig attribute*), 151  
 RELATIVE (*TiePieOscilloscopeTriggerLevelMode attribute*), 160  
 RELATIVE\_POSITION\_MOTOR (*ILS2T.ActionsPtp attribute*), 138  
 RELATIVE\_POSITION\_TARGET (*ILS2T.ActionsPtp attribute*), 138

```

remote() (Technix property), 149
remove_device() (DeviceSequenceMixin method), 70
require_block_measurement_support (TiePieDeviceConfig attribute), 154
required_keys() (AsyncCommunicationProtocolConfig class method), 12
required_keys() (CryLasAttenuatorConfig class method), 73
required_keys() (CryLasAttenuatorSerialCommunicationConfig class method), 74
required_keys() (CryLasLaserConfig class method), 78
required_keys() (CryLasLaserSerialCommunicationConfig class method), 80
required_keys() (EmptyConfig class method), 71, 166
required_keys() (FuGConfig class method), 87
required_keys() (FuGSerialCommunicationConfig class method), 95
required_keys() (HeinzingerConfig class method), 97
required_keys() (HeinzingerSerialCommunicationConfig class method), 101
required_keys() (ILS2TConfig class method), 141
required_keys() (ILS2TModbusTcpCommunicationConfig class method), 142
required_keys() (LJMCommunicationConfig class method), 15
required_keys() (LuminoxConfig class method), 145
required_keys() (LuminoxSerialCommunicationConfig class method), 148
required_keys() (MBW973Config class method), 109
required_keys() (MBW973SerialCommunicationConfig class method), 110
required_keys() (ModbusTcpCommunicationConfig class method), 17
required_keys() (NewportSMC100PPConfig class method), 119
required_keys() (NewportSMC100PPSerialCommunicationConfig class method), 123
required_keys() (OpcUaCommunicationConfig class method), 20
required_keys() (PfeifferTPGConfig class method), 128
required_keys() (PfeifferTPGSerialCommunicationConfig class method), 129
required_keys() (PSI9000Config class method), 84
required_keys() (PSI9000VisaCommunicationConfig class method), 85
required_keys() (RTO1024Config class method), 136
required_keys() (RTO1024VisaCommunicationConfig class method), 137
required_keys() (SerialCommunicationConfig class method), 23
required_keys() (SupercubeAOpcUaConfiguration class method), 50, 68
required_keys() (SupercubeBOpcUaConfiguration class method), 53
required_keys() (SupercubeConfiguration class method), 34, 57
required_keys() (SupercubeOpcUaCommunicationConfig class method), 34, 58
required_keys() (TechnixCommunicationConfig class method), 150
required_keys() (TechnixConfig class method), 151
required_keys() (TechnixSerialCommunicationConfig class method), 152
required_keys() (TechnixTelnetCommunicationConfig class method), 152
required_keys() (TelnetCommunicationConfig class method), 25
required_keys() (TiePieDeviceConfig class method), 154
required_keys() (VisaCommunicationConfig class method), 28
required_keys() (VisaDeviceConfig class method), 163
reset (BreakdownDetection attribute), 41, 60
RESET (FuGProbusIVCommands attribute), 90
reset () (FuGProbusIV method), 90
reset () (NewportSMC100PP method), 116
reset () (VisaDevice method), 162
reset_error () (ILS2T method), 140
reset_interface () (HeinzingerDI method), 98
resolution () (TiePieOscilloscopeConfig property), 159
response_sleep_time (CryLasAttenuatorConfig attribute), 73
RISING (TiePieOscilloscopeTriggerKind attribute), 160
RISING_OR_FALLING (TiePieOscilloscopeTriggerKind attribute), 160
RMS_CURRENT_LIMIT (NewportSMC100PP.MotorErrors attribute), 112
rpm_max_init (ILS2TConfig attribute), 142
rs485_address (NewportSMC100PPConfig attribute), 120
RTO1024 (class in hvl_ccb.dev.rs_rto1024), 130
RTO1024.TriggerModes (class in hvl_ccb.dev.rs_rto1024), 130
RTO1024Config (class in hvl_ccb.dev.rs_rto1024), 136
RTO1024Error, 136

```

RTO1024VisaCommunication (class *in* SerialCommunicationParity (class *in* hvl\_ccb.dev.rs\_rto1024), 136 hvl\_ccb.comm.serial), 23

RTO1024VisaCommunicationConfig (class *in* SerialCommunicationStopbits (class *in* hvl\_ccb.dev.rs\_rto1024), 136 hvl\_ccb.comm.serial), 23

run() (*ExperimentManager method*), 168

run\_continuous\_acquisition() (*RTO1024 method*), 133

run\_single\_acquisition() (*RTO1024 method*), 133

RUNNING (*ExperimentStatus attribute*), 168

## S

S (*LabJack.ThermocoupleType attribute*), 104

SA (*NewportConfigCommands attribute*), 111

safe\_ground\_enabled() (*TiePieOscilloscopeChannelConfig property*), 158

Safety (class *in* hvl\_ccb.dev.supercube.constants), 48

Safety (class *in* hvl\_ccb.dev.supercube2015.constants), 65

SafetyStatus (class *in* hvl\_ccb.dev.supercube.constants), 48

SafetyStatus (class *in* hvl\_ccb.dev.supercube2015.constants), 65

sample\_frequency() (*TiePieOscilloscopeConfig property*), 159

save\_configuration() (*RTO1024 method*), 133

save\_waveform\_history() (*RTO1024 method*), 133

SCALE (*ILS2TRegAddr attribute*), 143

ScalingFactorValueError, 143

screw\_scaling (*NewportSMC100PPConfig attribute*), 120

send\_command() (*NewportSMC100PPSerialCommunication method*), 122

send\_command() (*PfeifferTPGSerialCommunication method*), 128

send\_stop() (*NewportSMC100PPSerialCommunication method*), 122

Sensor\_error (*PfeifferTPG.SensorStatus attribute*), 125

Sensor\_off (*PfeifferTPG.SensorStatus attribute*), 125

sensor\_status (*LuminoxMeasurementType attribute*), 146

serial\_number (*LuminoxMeasurementType attribute*), 146

serial\_number (*TiePieDeviceConfig attribute*), 154

SerialCommunication (class *in* hvl\_ccb.comm.serial), 21

SerialCommunicationBytesize (class *in* hvl\_ccb.comm.serial), 22

SerialCommunicationConfig (class *in* hvl\_ccb.comm.serial), 22

SerialCommunicationIOError, 23

set\_acceleration() (*NewportSMC100PP method*), 116

set\_acquire\_length() (*RTO1024 method*), 133

set\_ain\_differential() (*LabJack method*), 105

set\_ain\_range() (*LabJack method*), 106

set\_ain\_resistance() (*LabJack method*), 106

set\_ain\_resolution() (*LabJack method*), 106

set\_ain\_thermocouple() (*LabJack method*), 106

set\_attenuation() (*CryLasAttenuator method*), 72

set\_cee16\_socket() (*Supercube2015Base method*), 55

set\_cee16\_socket() (*SupercubeBase method*), 31

set\_channel\_offset() (*RTO1024 method*), 133

set\_channel\_position() (*RTO1024 method*), 134

set\_channel\_range() (*RTO1024 method*), 134

set\_channel\_scale() (*RTO1024 method*), 134

set\_channel\_state() (*RTO1024 method*), 134

set\_current() (*HeinzingerDI method*), 98

set\_current() (*HeinzingerPNC method*), 100

set\_digital\_output() (*LabJack method*), 107

set\_display\_unit() (*PfeifferTPG method*), 126

set\_earthing\_manual() (*Supercube2015Base method*), 55

set\_full\_scale\_mbar() (*PfeifferTPG method*), 126

set\_full\_scale\_unitless() (*PfeifferTPG method*), 126

set\_init\_attenuation() (*CryLasAttenuator method*), 72

set\_init\_shutter\_status() (*CryLasLaser method*), 76

set\_jog\_speed() (*ILS2T method*), 140

set\_lower\_limits() (*PSI9000 method*), 82

set\_max\_acceleration() (*ILS2T method*), 140

set\_max\_deceleration() (*ILS2T method*), 140

set\_max\_rpm() (*ILS2T method*), 140

set\_measuring\_options() (*MBW973 method*), 108

set\_message\_board() (*SupercubeBase method*), 32

set\_motor\_configuration() (*NewportSMC100PP method*), 116

set\_negative\_software\_limit() (*NewportSMC100PP method*), 117

set\_number\_of\_recordings() (*HeinzingerDI method*), 99

set\_output() (*PSI9000 method*), 82

set\_positive\_software\_limit() (*NewportSMC100PP method*), 117

set\_pulse\_energy () (*CryLasLaser method*), 76  
 set\_ramp\_type () (*ILS2T method*), 140  
 set\_reference\_point () (*RTO1024 method*), 135  
 set\_register () (*FuGProbusV method*), 91  
 set\_remote\_control () (*Supercube2015Base method*), 56  
 set\_remote\_control () (*SupercubeBase method*), 32  
 set\_repetition\_rate () (*CryLasLaser method*), 76  
 set\_repetitions () (*RTO1024 method*), 135  
 set\_slope () (*Supercube2015WithFU method*), 67  
 set\_slope () (*SupercubeWithFU method*), 51  
 set\_status\_board () (*SupercubeBase method*), 32  
 set\_support\_output () (*Supercube2015Base method*), 56  
 set\_support\_output () (*SupercubeBase method*), 32  
 set\_support\_output\_impulse () (*Supercube2015Base method*), 56  
 set\_support\_output\_impulse () (*SupercubeBase method*), 32  
 set\_system\_lock () (*PSI9000 method*), 82  
 set\_t13\_socket () (*Supercube2015Base method*), 56  
 set\_t13\_socket () (*SupercubeBase method*), 33  
 set\_target\_voltage () (*Supercube2015WithFU method*), 67  
 set\_target\_voltage () (*SupercubeWithFU method*), 51  
 set\_transmission () (*CryLasAttenuator method*), 72  
 set\_trigger\_level () (*RTO1024 method*), 135  
 set\_trigger\_mode () (*RTO1024 method*), 135  
 set\_trigger\_source () (*RTO1024 method*), 135  
 set\_upper\_limits () (*PSI9000 method*), 82  
 set\_voltage () (*HeinzingerDI method*), 99  
 set\_voltage () (*HeinzingerPNC method*), 100  
 set\_voltage\_current () (*PSI9000 method*), 83  
 SETCURRENT (*FuGProbusVRegisterGroups attribute*), 93  
 setup (*Power attribute*), 47, 64  
 setvalue () (*FuGProbusVSetRegisters property*), 94  
 SETVOLTAGE (*FuGProbusVRegisterGroups attribute*), 93  
 SEVENBITS (*SerialCommunicationBytesize attribute*), 22  
 SHORT\_CIRCUIT (*NewportSMC100PP.MotorErrors attribute*), 112  
 SHUTDOWN\_CURRENT\_LIMIT (*PSI9000 attribute*), 81  
 SHUTDOWN\_VOLTAGE\_LIMIT (*PSI9000 attribute*), 81  
 ShutterStatus (*CryLasLaser attribute*), 75  
 ShutterStatus (*CryLasLaserConfig attribute*), 77  
 signal\_type () (*TiePieGeneratorConfig property*), 155  
 SINE (*TiePieGeneratorSignalType attribute*), 155  
 SingleCommDevice (*class in hvl\_ccb.dev.base*), 71  
 SIXBITS (*SerialCommunicationBytesize attribute*), 22  
 SIXTEEN (*HeinzingerConfig.RecordingsEnum attribute*), 96  
 SIXTEEN\_BIT (*TiePieOscilloscopeResolution attribute*), 160  
 SL (*NewportConfigCommands attribute*), 111  
 SN (*FuGReadbackChannels attribute*), 94  
 SOFTWARE\_INTERNAL\_SIXTY (*CryLasLaser.RepetitionRates attribute*), 75  
 SOFTWARE\_INTERNAL\_TEN (*CryLasLaser.RepetitionRates attribute*), 75  
 SOFTWARE\_INTERNAL\_TWENTY (*CryLasLaser.RepetitionRates attribute*), 75  
 software\_revision (*LuminoxMeasurementType attribute*), 146  
 SPACE (*SerialCommunicationParity attribute*), 23  
 SPECIALRAMPUPWARDS (*FuGRampModes attribute*), 94  
 spoll () (*VisaCommunication method*), 27  
 spoll\_handler () (*VisaDevice method*), 162  
 SQUARE (*TiePieGeneratorSignalType attribute*), 155  
 SR (*NewportConfigCommands attribute*), 111  
 srq\_mask () (*FuGProbusVConfigRegisters property*), 91  
 srq\_status () (*FuGProbusVConfigRegisters property*), 91  
 stage\_configuration (*NewportSMC100PPConfig attribute*), 120  
 start () (*CryLasAttenuator method*), 72  
 start () (*CryLasLaser method*), 76  
 start () (*Device method*), 69  
 start () (*DeviceSequenceMixin method*), 70  
 start () (*ExperimentManager method*), 168  
 start () (*FuG method*), 87  
 start () (*FuGProbusIV method*), 90  
 start () (*HeinzingerDI method*), 99  
 start () (*HeinzingerPNC method*), 100  
 start () (*ILS2T method*), 140  
 start () (*LabJack method*), 107  
 start () (*Luminox method*), 145  
 start () (*MBW973 method*), 108  
 start () (*NewportSMC100PP method*), 117  
 start () (*PfeifferTPG method*), 127  
 start () (*PSI9000 method*), 83  
 start () (*RTO1024 method*), 136  
 start () (*SingleCommDevice method*), 71  
 start () (*Supercube2015Base method*), 56  
 start () (*SupercubeBase method*), 33  
 start () (*Technix method*), 149  
 start () (*TiePieGeneratorMixin method*), 155  
 start () (*TiePieI2CHostMixin method*), 156

start () (*TiePieOscilloscope* method), 157  
 start () (*VisaDevice* method), 163  
 start\_control () (*MBW973* method), 108  
 start\_measurement () (*TiePieOscilloscope* method), 157  
 start\_polling () (*Poller* method), 161  
 STARTING (*ExperimentStatus* attribute), 168  
 States (*NewportSMC100PP* attribute), 113  
 status (*Safety* attribute), 48  
 status () (*EarthingStick* class method), 42  
 status () (*ExperimentManager* property), 168  
 status () (*FuGProbusVConfigRegisters* property), 91  
 status () (*FuGProbusVDORegisters* property), 92  
 status\_1 (*Door* attribute), 41  
 status\_1 (*EarthingRod* attribute), 41  
 status\_1 (*EarthingStick* attribute), 43  
 status\_1\_closed (*EarthingStick* attribute), 60  
 status\_1\_connected (*EarthingStick* attribute), 61  
 status\_1\_open (*EarthingStick* attribute), 61  
 status\_2 (*Door* attribute), 41  
 status\_2 (*EarthingRod* attribute), 41  
 status\_2 (*EarthingStick* attribute), 43  
 status\_2\_closed (*EarthingStick* attribute), 61  
 status\_2\_connected (*EarthingStick* attribute), 61  
 status\_2\_open (*EarthingStick* attribute), 61  
 status\_3 (*Door* attribute), 41  
 status\_3 (*EarthingRod* attribute), 41  
 status\_3 (*EarthingStick* attribute), 43  
 status\_3\_closed (*EarthingStick* attribute), 61  
 status\_3\_connected (*EarthingStick* attribute), 61  
 status\_3\_open (*EarthingStick* attribute), 61  
 status\_4 (*EarthingStick* attribute), 43  
 status\_4\_closed (*EarthingStick* attribute), 61  
 status\_4\_connected (*EarthingStick* attribute), 61  
 status\_4\_open (*EarthingStick* attribute), 61  
 status\_5 (*EarthingStick* attribute), 43  
 status\_5\_closed (*EarthingStick* attribute), 61  
 status\_5\_connected (*EarthingStick* attribute), 61  
 status\_5\_open (*EarthingStick* attribute), 61  
 status\_6 (*EarthingStick* attribute), 43  
 status\_6\_closed (*EarthingStick* attribute), 61  
 status\_6\_connected (*EarthingStick* attribute), 61  
 status\_6\_open (*EarthingStick* attribute), 61  
 status\_closed () (*EarthingStick* class method), 61  
 status\_connected () (*EarthingStick* class method), 61  
 status\_error (*Safety* attribute), 65  
 status\_green (*Safety* attribute), 65  
 status\_open () (*EarthingStick* class method), 61  
 status\_ready\_for\_red (*Safety* attribute), 65  
 status\_red (*Safety* attribute), 65  
 STATUSBYTE (*FuGReadbackChannels* attribute), 94  
 statuses () (*EarthingStick* class method), 43  
 stop (*Errors* attribute), 44, 62  
 stop () (*CryLasLaser* method), 77  
 stop () (*Device* method), 69  
 stop () (*DeviceSequenceMixin* method), 70  
 stop () (*ExperimentManager* method), 168  
 stop () (*FuGProbusIV* method), 90  
 stop () (*HeinzingerDI* method), 99  
 stop () (*ILS2T* method), 140  
 stop () (*LabJack* method), 107  
 stop () (*Luminox* method), 145  
 stop () (*MBW973* method), 108  
 stop () (*NewportSMC100PP* method), 117  
 stop () (*PfeifferTPG* method), 127  
 stop () (*PSI9000* method), 83  
 stop () (*RTO1024* method), 136  
 stop () (*SingleCommDevice* method), 71  
 stop () (*Supercube2015Base* method), 56  
 stop () (*SupercubeBase* method), 33  
 stop () (*Technix* method), 149  
 stop () (*TiePieGeneratorMixin* method), 155  
 stop () (*TiePieI2CHostMixin* method), 156  
 stop () (*TiePieOscilloscope* method), 157  
 stop () (*VisaDevice* method), 163  
 stop\_acquisition () (*RTO1024* method), 136  
 stop\_motion () (*NewportSMC100PP* method), 117  
 stop\_number (*Errors* attribute), 62  
 stop\_polling () (*Poller* method), 161  
 stopbits (*CryLasAttenuatorSerialCommunicationConfig* attribute), 74  
 stopbits (*CryLasLaserSerialCommunicationConfig* attribute), 80  
 stopbits (*FuGSerialCommunicationConfig* attribute), 95  
 stopbits (*HeinzingerSerialCommunicationConfig* attribute), 101  
 stopbits (*LuminoxSerialCommunicationConfig* attribute), 148  
 stopbits (*MBW973SerialCommunicationConfig* attribute), 110  
 stopbits (*NewportSMC100PPSerialCommunicationConfig* attribute), 123  
 stopbits (*PfeifferTPGSerialCommunicationConfig* attribute), 129  
 Stopbits (*SerialCommunicationConfig* attribute), 22  
 stopbits (*SerialCommunicationConfig* attribute), 23  
 streaming (*LuminoxOutputMode* attribute), 146  
 StrEnumBase (class in *hvl\_ccb.utils.enum*), 164  
 sub\_handler (*OpcUaCommunicationConfig* attribute), 20  
 sub\_handler (*SupercubeOpcUaCommunicationConfig* attribute), 34, 58  
 suitable\_range () (*TiePieOscilloscopeRange* static method), 159  
 Supercube2015Base (class in *hvl\_ccb.dev.supercube2015.base*), 54

Supercube2015WithFU	(class <i>hvl_ccb.dev.supercube2015.typ_a</i> ),	66	<i>in</i>	t13_2 ( <i>GeneralSockets attribute</i> ),	44, 62	
SupercubeAOpcUaCommunication	(class <i>hvl_ccb.dev.supercube.typ_a</i> ),	49	<i>in</i>	T13_SOCKET_PORTS	( <i>in module</i> <i>hvl_ccb.dev.supercube.constants</i> ),	48
SupercubeAOpcUaCommunication	(class <i>hvl_ccb.dev.supercube2015.typ_a</i> ),	67	<i>in</i>	T13_SOCKET_PORTS	( <i>in module</i> <i>hvl_ccb.dev.supercube2015.constants</i> ),	66
SupercubeAOpcUaConfiguration	(class <i>hvl_ccb.dev.supercube.typ_a</i> ),	49	<i>in</i>	T1MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeAOpcUaConfiguration	(class <i>hvl_ccb.dev.supercube2015.typ_a</i> ),	67	<i>in</i>	T200MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeB	(class <i>in hvl_ccb.dev.supercube.typ_b</i> ),	52	<i>in</i>	T20MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeBase	(class <i>hvl_ccb.dev.supercube.base</i> ),	29	<i>in</i>	T256US ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeBOpcUaCommunication	(class <i>hvl_ccb.dev.supercube.typ_b</i> ),	52	<i>in</i>	T4 ( <i>LabJack.DeviceType attribute</i> ),	103	
SupercubeBOpcUaConfiguration	(class <i>hvl_ccb.dev.supercube.typ_b</i> ),	52	<i>in</i>	T4 (LJMCommunicationConfig.DeviceType attribute),	14	
SupercubeBOpcUaConfiguration	(class <i>hvl_ccb.dev.supercube.base</i> ),	33	<i>in</i>	T40MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeConfiguration	(class <i>hvl_ccb.dev.supercube2015.base</i> ),	56	<i>in</i>	T4MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeEarthngStickOperationError,		34	<i>in</i>	T7 ( <i>LabJack.DeviceType attribute</i> ),	103	
SupercubeOpcEndpoint	(class <i>hvl_ccb.dev.supercube.constants</i> ),	48	<i>in</i>	T7 (LJMCommunicationConfig.DeviceType attribute),	14	
SupercubeOpcEndpoint	(class <i>hvl_ccb.dev.supercube2015.constants</i> ),	66	<i>in</i>	T7_PRO ( <i>LabJack.DeviceType attribute</i> ),	103	
SupercubeOpcUaCommunication	(class <i>hvl_ccb.dev.supercube.base</i> ),	34	<i>in</i>	T7_PRO (LJMCommunicationConfig.DeviceType attribute),	15	
SupercubeOpcUaCommunication	(class <i>hvl_ccb.dev.supercube2015.base</i> ),	57	<i>in</i>	T800MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeOpcUaCommunicationConfig	(class <i>in hvl_ccb.dev.supercube.base</i> ),	34	<i>in</i>	T80MS ( <i>FuGMonitorModes attribute</i> ),	89	
SupercubeOpcUaCommunicationConfig	(class <i>in hvl_ccb.dev.supercube2015.base</i> ),	57	<i>in</i>	target_pulse_energy () ( <i>CryLasLaser property</i> ),	77	
SupercubeSubscriptionHandler	(class <i>hvl_ccb.dev.supercube.base</i> ),	34	<i>in</i>	TCP (LJMCommunicationConfig.ConnectionType attribute),	14	
SupercubeSubscriptionHandler	(class <i>hvl_ccb.dev.supercube2015.base</i> ),	58	<i>in</i>	TCPIP_INSTR (VisaCommunicationConfig.InterfaceType attribute),	27	
SupercubeWithFU	(class <i>hvl_ccb.dev.supercube.typ_a</i> ),	50	<i>in</i>	TCPIP_SOCKET (VisaCommunicationConfig.InterfaceType attribute),	27	
switch_to_operate ( <i>Safety attribute</i> ),	48			TEC1 ( <i>CryLasLaser.AnswersStatus attribute</i> ),	75	
switch_to_ready ( <i>Safety attribute</i> ),	48			TEC2 ( <i>CryLasLaser.AnswersStatus attribute</i> ),	75	
switchto_green ( <i>Safety attribute</i> ),	65			Technix (class <i>in hvl_ccb.dev.technix</i> ),	148	
switchto_operate ( <i>Safety attribute</i> ),	65			TechnixCommunication	(class <i>in hvl_ccb.dev.technix</i> ),	149
switchto_ready ( <i>Safety attribute</i> ),	65			TechnixCommunicationConfig	(class <i>in hvl_ccb.dev.technix</i> ),	149
SyncCommunicationProtocol	(class <i>hvl_ccb.comm.base</i> ),	12	<i>in</i>	TechnixConfig (class <i>in hvl_ccb.dev.technix</i> ),	150	
SyncCommunicationProtocolConfig	(class <i>in</i> <i>hvl_ccb.comm.base</i> ),	12	<i>in</i>	TechnixError,	151	
<b>T</b>				TechnixSerialCommunication	(class <i>in hvl_ccb.dev.technix</i> ),	151
T ( <i>LabJack.ThermocoupleType attribute</i> ),	104			TechnixSerialCommunicationConfig	(class <i>in hvl_ccb.dev.technix</i> ),	151
t13_1 ( <i>GeneralSockets attribute</i> ),	44, 62			TechnixStatusByte	(class <i>in hvl_ccb.dev.technix</i> ),	152
				TechnixTelnetCommunication	(class <i>in hvl_ccb.dev.technix</i> ),	152
				TechnixTelnetCommunicationConfig	(class <i>in hvl_ccb.dev.technix</i> ),	152
				TelnetCommunication	(class <i>in hvl_ccb.comm.telnet</i> ),	24
				TelnetCommunicationConfig	(class <i>in hvl_ccb.comm.telnet</i> ),	25
				TelnetError,	25	
				TEMP (ILS2TRegAddr attribute),	143	

temperature\_sensor (*LuminoxMeasurementType attribute*), 146  
TEN (*LabJack.AInRange attribute*), 103  
TEN (*LabJack.CalMicroAmpere attribute*), 103  
terminator (*AsyncCommunicationProtocolConfig attribute*), 12  
terminator (*CryLasAttenuatorSerialCommunicationConfig attribute*), 74  
terminator (*CryLasLaserSerialCommunicationConfig attribute*), 80  
TERMINATOR (*FuGProbusIVCommands attribute*), 90  
terminator (*FuGSerialCommunicationConfig attribute*), 95  
terminator (*HeinzingerSerialCommunicationConfig attribute*), 101  
terminator (*LuminoxSerialCommunicationConfig attribute*), 148  
terminator (*MBW973SerialCommunicationConfig attribute*), 110  
terminator (*NewportSMC100PPSerialCommunicationConfig attribute*), 123  
terminator (*PfeifferTPGSerialCommunicationConfig attribute*), 129  
terminator (*TechnixCommunicationConfig attribute*), 150  
terminator() (*FuGProbusVConfigRegisters property*), 91  
terminator\_str() (*SerialCommunicationConfig method*), 23  
TiePieDeviceConfig (*class in hvl\_ccb.dev.tiepie*), 153  
TiePieDeviceType (*class in hvl\_ccb.dev.tiepie*), 154  
TiePieError, 154  
TiePieGeneratorConfig (*class in hvl\_ccb.dev.tiepie*), 154  
TiePieGeneratorMixin (*class in hvl\_ccb.dev.tiepie*), 155  
TiePieGeneratorSignalType (*class in hvl\_ccb.dev.tiepie*), 155  
TiePieHS5 (*class in hvl\_ccb.dev.tiepie*), 155  
TiePieHS6 (*class in hvl\_ccb.dev.tiepie*), 156  
TiePieI2CHostConfig (*class in hvl\_ccb.dev.tiepie*), 156  
TiePieI2CHostMixin (*class in hvl\_ccb.dev.tiepie*), 156  
TiePieOscilloscope (*class in hvl\_ccb.dev.tiepie*), 156  
TiePieOscilloscopeAutoResolutionModes (*class in hvl\_ccb.dev.tiepie*), 157  
TiePieOscilloscopeChannelConfig (*class in hvl\_ccb.dev.tiepie*), 158  
TiePieOscilloscopeChannelCoupling (*class in hvl\_ccb.dev.tiepie*), 158  
TiePieOscilloscopeConfig (*class in hvl\_ccb.dev.tiepie*), 159  
TiePieOscilloscopeRange (*class in hvl\_ccb.dev.tiepie*), 159  
TiePieOscilloscopeResolution (*class in hvl\_ccb.dev.tiepie*), 159  
TiePieOscilloscopeTriggerKind (*class in hvl\_ccb.dev.tiepie*), 160  
TiePieOscilloscopeTriggerLevelMode (*class in hvl\_ccb.dev.tiepie*), 160  
TiePieWS5 (*class in hvl\_ccb.dev.tiepie*), 160  
timeout (*CryLasAttenuatorSerialCommunicationConfig attribute*), 74  
timeout (*CryLasLaserSerialCommunicationConfig attribute*), 80  
timeout (*FuGSerialCommunicationConfig attribute*), 95  
timeout (*HeinzingerSerialCommunicationConfig attribute*), 102  
timeout (*LuminoxSerialCommunicationConfig attribute*), 148  
timeout (*MBW973SerialCommunicationConfig attribute*), 110  
timeout (*NewportSMC100PPSerialCommunicationConfig attribute*), 123  
timeout (*PfeifferTPGSerialCommunicationConfig attribute*), 129  
timeout (*SerialCommunicationConfig attribute*), 23  
timeout (*TelnetCommunicationConfig attribute*), 25  
timeout (*VisaCommunicationConfig attribute*), 28  
Torr (*PfeifferTPG.PressureUnits attribute*), 124  
TPG25xA (*PfeifferTPGConfig.Model attribute*), 127  
TPGx6x (*PfeifferTPGConfig.Model attribute*), 127  
TPR (*PfeifferTPG.SensorTypes attribute*), 125  
transmission() (*CryLasAttenuator property*), 72  
TRIANGLE (*TiePieGeneratorSignalType attribute*), 155  
trigger\_enabled() (*TiePieOscilloscopeChannelConfig property*), 158  
trigger\_hysteresis() (*TiePieOscilloscopeChannelConfig property*), 158  
trigger\_kind() (*TiePieOscilloscopeChannelConfig property*), 158  
trigger\_level() (*TiePieOscilloscopeChannelConfig property*), 158  
trigger\_level\_mode() (*TiePieOscilloscopeChannelConfig property*), 158  
trigger\_timeout() (*TiePieOscilloscopeConfig property*), 159  
triggered (*BreakdownDetection attribute*), 41, 60  
TWELVE\_BIT (*TiePieOscilloscopeResolution attribute*), 160  
TWENTY\_VOLT (*TiePieOscilloscopeRange attribute*), 159  
TWO (*HeinzingerConfig.RecordingsEnum attribute*), 96  
TWO (*SerialCommunicationStopbits attribute*), 24

TWO\_HUNDRED (*LabJack.CalMicroAmpere attribute*), 103  
 TWO\_HUNDRED\_MILLI\_VOLT (*TiePieOscilloscopeRange attribute*), 159  
 TWO\_VOLT (*TiePieOscilloscopeRange attribute*), 159

**U**

Underrange (*PfeifferTPG.SensorStatus attribute*), 125  
 unit (*ILS2TModbusTcpCommunicationConfig attribute*), 142  
 unit (*ModbusTcpCommunicationConfig attribute*), 17  
 unit () (*PfeifferTPG property*), 127  
 unit\_current () (*HeinzingerPNC property*), 100  
 unit\_voltage () (*HeinzingerPNC property*), 100  
 UNKNOWN (*HeinzingerDI.OutputStatus attribute*), 97  
 UNKNOWN (*HeinzingerPNC.UnitCurrent attribute*), 99  
 UNKNOWN (*HeinzingerPNC.UnitVoltage attribute*), 99  
 UNKNOWN (*TiePieGeneratorSignalType attribute*), 155  
 UNKNOWN (*TiePieOscilloscopeAutoResolutionModes attribute*), 158  
 UNKNOWN (*TiePieOscilloscopeTriggerLevelMode attribute*), 160  
 UNREADY\_INACTIVE (*CryLasLaser.LaserStatus attribute*), 75  
 update\_laser\_status () (*CryLasLaser method*), 77  
 update\_period (*OpcUaCommunicationConfig attribute*), 20  
 update\_repetition\_rate () (*CryLasLaser method*), 77  
 update\_shutter\_status () (*CryLasLaser method*), 77  
 update\_target\_pulse\_energy () (*CryLasLaser method*), 77  
 UpdateEspStageInfo (*New-portSMC100PPConfig.EspStageConfig attribute*), 118  
 USB (*LJMCommunicationConfig.ConnectionType attribute*), 14  
 user\_position\_offset (*New-portSMC100PPConfig attribute*), 120  
 user\_steps () (*ILS2T method*), 141

**V**

V (*HeinzingerPNC.UnitVoltage attribute*), 99  
 VA (*NewportConfigCommands attribute*), 111  
 value () (*FuGProbusVMonitorRegisters property*), 92  
 value () (*LabJack.AInRange property*), 103  
 value\_raw () (*FuGProbusVMonitorRegisters property*), 93  
 ValueEnum (*class in hvl\_ccb.utils.enum*), 164  
 VB (*NewportConfigCommands attribute*), 111  
 velocity (*NewportSMC100PPConfig attribute*), 120

visa\_backend (*VisaCommunicationConfig attribute*), 28  
 VisaCommunication (*class in hvl\_ccb.comm.visa*), 26  
 VisaCommunicationConfig (*class in hvl\_ccb.comm.visa*), 27  
 VisaCommunicationConfig.InterfaceType (*class in hvl\_ccb.comm.visa*), 27  
 VisaCommunicationError, 28  
 VisaDevice (*class in hvl\_ccb.dev.visa*), 162  
 VisaDeviceConfig (*class in hvl\_ccb.dev.visa*), 163  
 VOLT (*ILS2TRegAddr attribute*), 143  
 Volt (*PfeifferTPG.PressureUnits attribute*), 124  
 VOLTAGE (*FuGProbusIVCommands attribute*), 90  
 VOLTAGE (*FuGReadbackChannels attribute*), 94  
 voltage () (*FuG property*), 87  
 voltage () (*Technix property*), 149  
 voltage\_lower\_limit (*PSI9000Config attribute*), 84  
 voltage\_max (*Power attribute*), 47, 64  
 voltage\_monitor () (*FuG property*), 87  
 voltage\_primary (*Power attribute*), 47, 64  
 voltage\_regulation () (*Technix property*), 149  
 voltage\_slope (*Power attribute*), 47, 64  
 voltage\_target (*Power attribute*), 47, 64  
 voltage\_upper\_limit (*PSI9000Config attribute*), 84

**W**

WAIT\_AFTER\_WRITE (*VisaCommunication attribute*), 26  
 wait\_for\_polling\_result () (*Poller method*), 161  
 wait\_operation\_complete () (*VisaDevice method*), 163  
 wait\_sec\_initialisation (*PSI9000Config attribute*), 84  
 wait\_sec\_max\_disable (*ILS2TConfig attribute*), 142  
 wait\_sec\_post\_absolute\_position (*ILS2TConfig attribute*), 142  
 wait\_sec\_post\_activate (*LuminoxConfig attribute*), 145  
 wait\_sec\_post\_CANNOT\_DISABLE (*ILS2TConfig attribute*), 142  
 wait\_sec\_post\_enable (*ILS2TConfig attribute*), 142  
 wait\_sec\_post\_relative\_step (*ILS2TConfig attribute*), 142  
 wait\_sec\_read\_text\_nonempty (*AsyncCommunicationProtocolConfig attribute*), 12  
 wait\_sec\_read\_text\_nonempty (*FuGSerial-CommunicationConfig attribute*), 96

**Z**  
 wait\_sec\_read\_text\_nonempty (*HeinzingerSerialCommunicationConfig attribute*), 102  
 wait\_sec\_retry\_get\_device (*TiePieDeviceConfig attribute*), 154  
 wait\_sec\_settings\_effect (*PSI9000Config attribute*), 84  
 wait\_sec\_stop\_commands (*FuGConfig attribute*), 88  
 wait\_sec\_stop\_commands (*HeinzingerConfig attribute*), 97  
 wait\_sec\_system\_lock (*PSI9000Config attribute*), 84  
 wait\_sec\_trials\_activate (*LuminoxConfig attribute*), 145  
 wait\_timeout\_retry\_sec (*OpcUaCommunicationConfig attribute*), 20  
 wait\_until\_motor\_initialized() (*NewportSMC100PP method*), 118  
 wait\_until\_ready() (*CryLasLaser method*), 77  
 warning (*Errors attribute*), 44  
 WIFI (*LJMCommunicationConfig.ConnectionType attribute*), 14  
 wrap\_libtiepie\_exception() (*in module hvl\_ccb.dev.tiepie*), 161  
 write() (*AsyncCommunicationProtocol method*), 10  
 write() (*MBW973 method*), 108  
 write() (*OpcUaCommunication method*), 19  
 write() (*Supercube2015Base method*), 56  
 write() (*SupercubeBase method*), 33  
 write() (*VisaCommunication method*), 27  
 write\_absolute\_position() (*ILS2T method*), 141  
 write\_bytes() (*AsyncCommunicationProtocol method*), 11  
 write\_bytes() (*SerialCommunication method*), 21  
 write\_bytes() (*TelnetCommunication method*), 25  
 write\_name() (*LJMCommunication method*), 14  
 write\_names() (*LJMCommunication method*), 14  
 write\_registers() (*ModbusTcpCommunication method*), 16  
 write\_relative\_step() (*ILS2T method*), 141  
 write\_termination (*VisaCommunicationConfig attribute*), 28  
 write\_text() (*AsyncCommunicationProtocol method*), 11  
 WRONG\_ESP\_STAGE (*NewportSMC100PP.MotorErrors attribute*), 112

**X**

x\_stat() (*FuGProbusVDIRegisters property*), 92  
 XOUTPUTS (*FuGProbusIVCommands attribute*), 91

**Y**

YES (*FuGDigitalVal attribute*), 88