
HVL Common Code Base Documentation

Release 0.8.3

Mikolaj Rybiński, David Graber, Henrik Menne, Alise Chachereau,

Sep 27, 2021

CONTENTS:

1	HVL Common Code Base	1
1.1	Features	1
1.2	Documentation	5
1.3	Credits	5
2	Installation	7
2.1	Stable release	7
2.2	From sources	7
2.3	Additional system libraries	8
3	Usage	9
4	API Documentation	11
4.1	hvl_ccb	11
5	Contributing	187
5.1	Types of Contributions	187
5.2	Get Started!	188
5.3	Merge Request Guidelines	189
5.4	Tips	189
5.5	Deploying	190
6	Credits	191
6.1	Maintainers	191
6.2	Authors	191
6.3	Contributors	191
7	History	193
7.1	0.8.3 (2021-09-27)	193
7.2	0.8.2 (2021-08-27)	193
7.3	0.8.1 (2021-08-13)	193
7.4	0.8.0 (2021-07-02)	193
7.5	0.7.1 (2021-06-04)	194
7.6	0.7.0 (2021-05-25)	194
7.7	0.6.1 (2021-05-08)	194
7.8	0.6.0 (2021-04-23)	195
7.9	0.5.0 (2020-11-11)	195
7.10	0.4.0 (2020-07-16)	196
7.11	0.3.5 (2020-02-18)	196
7.12	0.3.4 (2019-12-20)	197
7.13	0.3.3 (2019-05-08)	197

7.14	0.3.2 (2019-05-08)	197
7.15	0.3.1 (2019-05-02)	197
7.16	0.3 (2019-05-02)	197
7.17	0.2.1 (2019-04-01)	198
7.18	0.2.0 (2019-03-31)	198
7.19	0.1.0 (2019-02-06)	198
8	Indices and tables	199
	Python Module Index	201
	Index	203

**CHAPTER
ONE**

HVL COMMON CODE BASE

Python common code base to control devices high voltage research devices, in particular, as used in Christian Franck's High Voltage Lab (HVL), D-ITET, ETH.

- Free software: GNU General Public License v3
- Copyright (c) 2019-2021 ETH Zurich, SIS ID and HVL D-ITET

1.1 Features

For managing multi-device experiments instantiate the `ExperimentManager` utility class.

1.1.1 Devices

The devices wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, various settings and options enumerations, as well as start/stop methods. Currently, wrappers to control the following devices are available:

Function/Type	Devices
Data acquisition	LabJack (T4, T7, T7-PRO; requires LJM Library) Pico Technology PT-104 Platinum Resistance Data Logger (requires PicoSDK/libusbpt104)
Digital IO	LabJack (T4, T7, T7-PRO; requires LJM Library)
Experiment control	HVL Supercube with and without Frequency Converter
Gas Analyser	MBW 973-SF6 gas dew point mirror analyzer Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges SST Luminox oxygen sensor
I2C host	TiePie (HS5, WS5; requires LibTiePie SDK)
Laser	CryLaS pulsed laser CryLaS laser attenuator
Oscilloscope	Rhode & Schwarz RTO 1024 TiePie (HS5, HS6, WS5; requires LibTiePie SDK)
Power supply	Elektro-Automatik PSI9000 FuG Elektronik Heinzinger PNC Technix capacitor charger
Stepper motor drive	Newport SMC100PP Schneider Electric ILS2T
Temperature control	Lauda PRO RP 245 E circulation thermostat
Waveform generator	TiePie (HS5, WS5; requires LibTiePie SDK)

Each device uses at least one standardised communication protocol wrapper.

1.1.2 Communication protocols

In `hvl_ccb` by “communication protocol” we mean different levels of communication standards, from the low level actual communication protocols like serial communication to application level interfaces like VISA TCP standard. There are also devices in `hvl_ccb` that use dummy communication protocol concept; this is because these devices build on proprietary vendor libraries that communicate with vendor devices, like in case of the TiePie devices.

The communication protocol wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, as well as open/close, and read/write/query methods. Currently, wrappers to use the following communication protocols are available:

Communication protocol	Devices using
Modbus TCP	Schneider Electric ILS2T stepper motor drive
OPC UA	HVL Supercube with and without Frequency Converter
Serial	CryLaS pulsed laser and laser attenuator FuG Elektronik power supply (e.g. capacitor charger HCK) using the Probus V protocol Heinzinger PNC power supply using Heinzinger Digital Interface I/II SST Luminox oxygen sensor MBW 973-SF6 gas dew point mirror analyzer Newport SMC100PP single axis driver for 2-phase stepper motors Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges Technix capacitor charger
TCP	Lauda PRO RP 245 E circulation thermostat
Telnet	Technix capacitor charger
VISA TCP	Elektro-Automatik PSI9000 DC power supply Rhode & Schwarz RTO 1024 oscilloscope
<i>proprietary</i>	LabJack (T4, T7, T7-PRO) devices, which communicate via LJM Library Pico Technology PT-104 Platinum Resistance Data Logger, which communicate via PicoSDK/libusbpt104 TiePie (HS5, HS6, WS5) oscilloscopes, generators and I2C hosts, which communicate via LibTiePie SDK

1.2 Documentation

Note: if you're planning to contribute to the hvl_ccb project do read beforehand the **Contributing** section in the HVL CCB documentation.

Do either:

- read [HVL CCB documentation at RTD](#),

or

- build and read HVL CCB documentation locally; install first [Graphviz](#) (make sure to have the `dot` command in the executable search path) and the Python build requirements for documentation:

```
$ pip install docs/requirements.txt
```

and then either on Windows in Git BASH run:

```
$ ./make.sh docs
```

or from any other shell with GNU Make installed run:

```
$ make docs
```

The target index HTML ("docs/_build/html/index.html") should open automatically in your Web browser.

1.3 Credits

This package was created with [Cookiecutter](#) and the [audreyr/cookiecutter-pypackage](#) project template.

INSTALLATION

2.1 Stable release

To install HVL Common Code Base, run this command in your terminal:

```
$ pip install hvl_ccb
```

To install HVL Common Code Base with optional Python libraries that require manual installations of additional system libraries, you need to specify on installation extra requirements corresponding to these controllers. For instance, to install Python requirements for LabJack and TiePie devices, run:

```
$ pip install "hvl_ccb[tiepie,labjack]"
```

See below for the info about additional system libraries and the corresponding extra requirements.

To install all extra requirements run:

```
$ pip install "hvl_ccb[all]"
```

This is the preferred method to install HVL Common Code Base, as it will always install the most recent stable release.

If you don't have `pip` installed, this [Python installation guide](#) can guide you through the process.

2.2 From sources

The sources for HVL Common Code Base can be downloaded from the [GitLab](#) repo.

You can either clone the repository:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

Or download the [tarball](#):

```
$ curl -OL https://gitlab.com/ethz_hvl/hvl_ccb/-/archive/master/hvl_ccb.tar.gz
```

Once you have a copy of the source, you can install it with:

```
$ pip install .
```

2.3 Additional system libraries

If you have installed *hvl_ccb* with any of the extra features corresponding to device controllers, you must additionally install respective system library; these are:

Extra feature	Additional system library
labjack	LJM Library
picotech	PicoSDK (Windows) / libusbpt104 (Ubuntu/Debian)
tiepie	LibTiePie SDK

For more details on installation of the libraries see docstrings of the corresponding `hvl_ccb` modules.

**CHAPTER
THREE**

USAGE

To use HVL Common Code Base in a project:

```
import hvl_ccb
```


API DOCUMENTATION


4.1 hvl_ccb

4.1.1 Subpackages

[hvl_ccb.comm](#)

Submodules

[hvl_ccb.comm.base](#)

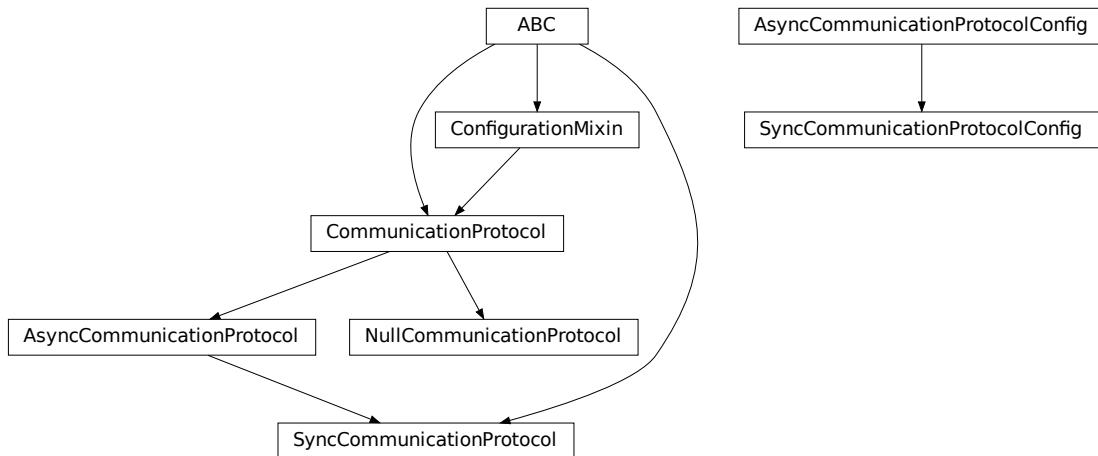

Module with base classes for communication protocols.

class AsyncCommunicationProtocol(config)
Bases: [hvl_ccb.comm.base.CommunicationProtocol](#)

Abstract base class for asynchronous communication protocols

static config_cls() → Type[[hvl_ccb.comm.base.AsyncCommunicationProtocolConfig](#)]
Return the default configdataclass class.

Returns a reference to the default configdataclass class

read() → str

Read a single line of text as *str* from the communication.

Returns text as *str* including the terminator, which can also be empty “”

read_all(*n_attempts_max*: Optional[int] = None, *attempt_interval_sec*: Optional[Union[int, float]] = None) → Optional[str]

Read all lines of text from the connection till nothing is left to read.

Parameters

- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt_interval_sec** – time between the reading attempts

Returns A multi-line *str* including the terminator internally

abstract read_bytes() → bytes

Read a single line as *bytes* from the communication.

This method uses *self.access_lock* to ensure thread-safety.

Returns a single line as *bytes* containing the terminator, which can also be empty b””

read_nonempty(*n_attempts_max*: Optional[int] = None, *attempt_interval_sec*: Optional[Union[int, float]] = None) → Optional[str]

Try to read a non-empty single line of text as *str* from the communication. If the host does not reply or reply with white space only, it will return None.

Returns a non-empty text as a *str* or *None* in case of an empty string

Parameters

- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt_interval_sec** – time between the reading attempts

read_text() → str

Read one line of text from the serial port. The input buffer may hold additional data afterwards, since only one line is read.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Returns String read from the serial port; “” if there was nothing to read.

Raises *SerialCommunicationIOError* – when communication port is not opened

read_text_nonempty(*n_attempts_max*: Optional[int] = None, *attempt_interval_sec*: Optional[Union[int, float]] = None) → Optional[str]

Reads from the serial port, until a non-empty line is found, or the number of attempts is exceeded.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Attention: in contrast to *read_text*, the returned answer will be stripped of a whitespace newline terminator at the end, if such terminator is set in the initial configuration (default).

Parameters

- **n_attempts_max** – maximum number of read attempts
- **attempt_interval_sec** – time between the reading attempts

Returns String read from the serial port; “” if number of attempts is exceeded or serial port is not opened.

write(*text*: str)

Write text as *str* to the communication.

Parameters `text` – test as a `str` to be written

abstract `write_bytes(data: bytes) → int`
 Write data as `bytes` to the communication.
 This method uses `self.access_lock` to ensure thread-safety.

Parameters `data` – data as `bytes`-string to be written

Returns number of bytes written

`write_text(text: str)`
 Write text to the serial port. The text is encoded and terminated by the configured terminator.
 NOTE: backward-compatibility proxy for `read` method; to be removed in v1.0

Parameters `text` – Text to send to the port.

Raises `SerialCommunicationIOError` – when communication port is not opened

```
class AsyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10)
```

Bases: `object`

Base configuration data class for asynchronous communication protocols

`clean_values()`

`default_n_attempts_read_text_nonempty: int = 10`
 default number of attempts to read a non-empty text

`encoding: str = 'utf-8'`
 Standard encoding of the connection. Typically this is `utf-8`, but can also be `latin-1` or something from here: <https://docs.python.org/3/library/codecs.html#standard-encodings>

`encoding_error_handling: str = 'strict'`
 Encoding error handling scheme as defined here: <https://docs.python.org/3/library/codecs.html#error-handlers> By default strict error handling that raises `UnicodeError`.

`force_value(fieldname, value)`
 Forces a value to a dataclass field despite the class being frozen.
 NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- `fieldname` – name of the field
- `value` – value to assign

`is_configdataclass = True`

`classmethod keys() → Sequence[str]`
 Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

`classmethod optional_defaults() → Dict[str, object]`
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

```
classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

terminator: bytes = b'\r\n'
    The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations.
    This defines the end of a single line.

wait_sec_read_text_nonempty: Union[int, float] = 0.5
    time to wait between attempts of reading a non-empty text

class CommunicationProtocol(config)
    Bases: hvl\_ccb.configuration.ConfigurationMixin, abc.ABC
    Communication protocol abstract base class.

    Specifies the methods to implement for communication protocol, as well as implements some default settings and checks.

access_lock
    Access lock to use with context manager when accessing the communication protocol (thread safety)

abstract close()
    Close the communication protocol

abstract open()
    Open communication protocol

class NullCommunicationProtocol(config)
    Bases: hvl\_ccb.comm.base.CommunicationProtocol

    Communication protocol that does nothing.

close() → None
    Void close function.

static config_cls() → Type[hvl\_ccb.configuration.EmptyConfig]
    Empty configuration

    Returns EmptyConfig

open() → None
    Void open function.

class SyncCommunicationProtocol(config)
    Bases: hvl\_ccb.comm.base.AsyncCommunicationProtocol, abc.ABC
    Abstract base class for synchronous communication protocols with query()

static config_cls() → Type[hvl\_ccb.comm.base.SyncCommunicationProtocolConfig]
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class


query(command: str) → Optional[str]
    Send a command to the interface and handle the status message. Eventually raises an exception.

    Parameters command – Command to send

    Returns Answer from the interface, which can be None instead of an empty reply
```

```
class SyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                      encoding_error_handling: str = 'strict',
                                      wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                      default_n_attempts_read_text_nonempty: int = 10)
Bases: hvl_ccb.comm.base.AsyncCommunicationProtocolConfig
```

hvl_ccb.comm.labjack_ljm

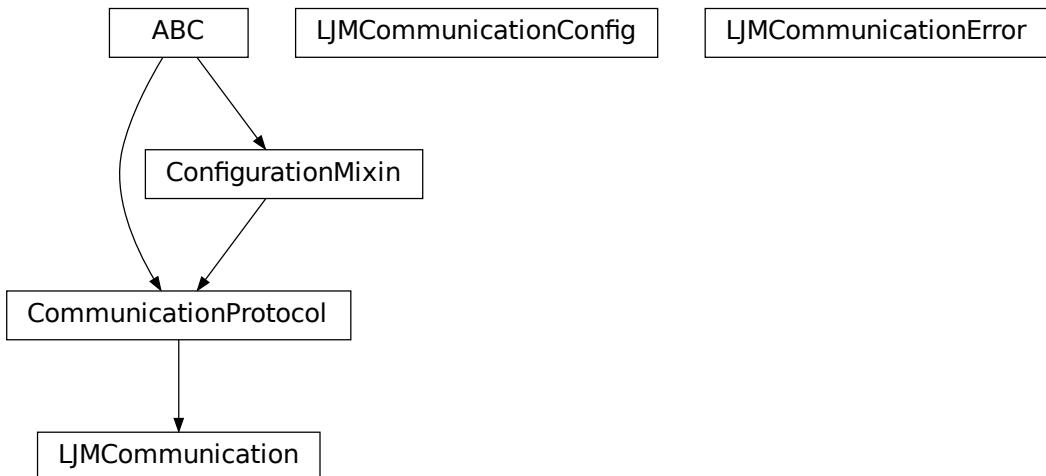

Communication protocol for LabJack using the LJM Library. Originally developed and tested for LabJack T7-PRO. Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

```
class LJMCommunication(configuration)
Bases: hvl_ccb.comm.base.CommunicationProtocol

Communication protocol implementing the LabJack LJM Library Python wrapper.

close() → None
    Close the communication port.

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

property is_open: bool
    Flag indicating if the communication port is open.

    Returns True if the port is open, otherwise False

open() → None
    Open the communication port.
```

read_name(*names: str, return_num_type: Type[numbers.Real] = <class 'float'>) → Union[numbers.Real, Sequence[numbers.Real]]

Read one or more input numeric values by name.

Parameters

- **names** – one or more names to read out from the LabJack
- **return_num_type** – optional numeric type specification for return values; by default *float*.

Returns answer of the LabJack, either single number or multiple numbers in a sequence, respectively, when one or multiple names to read were given

Raises **TypeError** – if read value of type not compatible with *return_num_type*

write_name(name: str, value: numbers.Real) → None

Write one value to a named output.

Parameters

- **name** – String or with name of LabJack IO
- **value** – is the value to write to the named IO port

write_names(name_value_dict: Dict[str, numbers.Real]) → None

Write more than one value at once to named outputs.

Parameters **name_value_dict** – is a dictionary with string names of LabJack IO as keys and corresponding numeric values

```
class LJMCommunicationConfig(device_type: Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY',
                               connection_type: Union[str,
                               hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType] =
                               'ANY', identifier: str = 'ANY')
```

Bases: object

Configuration dataclass for [LJMCommunication](#).

```
class ConnectionType(value=<no_arg>, names=None, module=None, type=None, start=1,
                      boundary=None)
```

Bases: [hvl_ccb.utils.enum.AutoNumberNameEnum](#)

LabJack connection type.

ANY = 1

ETHERNET = 4

TCP = 3

USB = 2

WIFI = 5

```
class DeviceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: [hvl_ccb.utils.enum.AutoNumberNameEnum](#)

LabJack device types.

Can be also looked up by ambiguous Product ID (*p_id*) or by instance name: `python LabJackDeviceType(4)` is `LabJackDeviceType('T4')`

ANY = 1

T4 = 2

T7 = 3

```
T7_PRO = 4
```

classmethod `get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
List[hvl_ccb._dev.labjack.DeviceType]]`

Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

Parameters `p_id` – Product ID of a LabJack device

Returns Instance or list of instances of `LabJackDeviceType`

Raises `ValueError` – when Product ID is unknown

clean_values() → None

Performs value checks on `device_type` and `connection_type`.

connection_type: `Union[str,
hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType] = 'ANY'`

Can be either string or of enum `ConnectionType`.

device_type: `Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY'`

Can be either string ‘ANY’, ‘T7_PRO’, ‘T7’, ‘T4’, or of enum `DeviceType`.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- `fieldname` – name of the field
- `value` – value to assign

identifier: str = 'ANY'

The identifier specifies information for the connection to be used. This can be an IP address, serial number, or device name. See the LabJack docs (<https://labjack.com/support/software/api/ljm/function-reference/ljmopens/identifier-parameter>) for more information.

is_configdataclass = True

classmethod `keys() → Sequence[str]`

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod `optional_defaults() → Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod `required_keys() → Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns a list of strings containing all required keys.

exception LJMCommunicationError

Bases: `Exception`

Errors coming from LJMCommunication.

hvl_ccb.comm.modbus_tcp

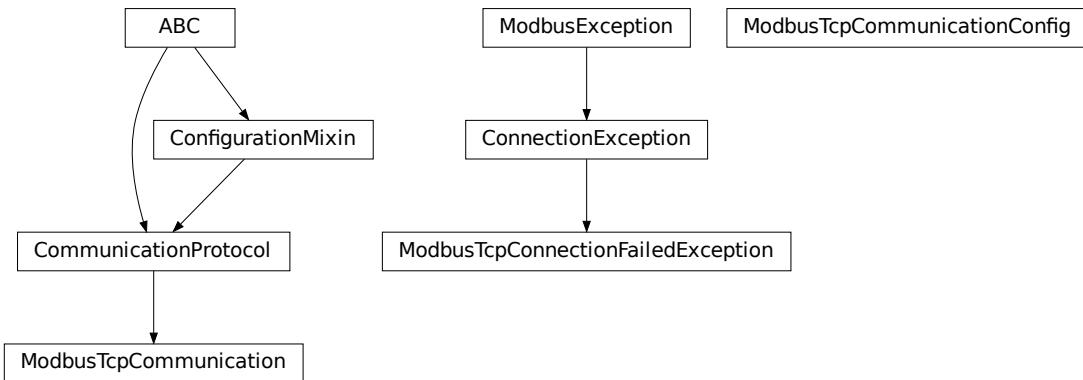

Communication protocol for modbus TCP ports. Makes use of the `pymodbus` library.

class ModbusTcpCommunication(configuration)
Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Implements the Communication Protocol for modbus TCP.

close()

Close the Modbus TCP connection.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

open() → None

Open the Modbus TCP connection.

Raises `ModbusTcpConnectionFailedException` – if the connection fails.

read_holding_registers(address: int, count: int) → List[int]

Read specified number of register starting with given address and return the values from each register.

Parameters

- **address** – address of the first register
- **count** – count of registers to read

Returns list of `int` values

read_input_registers(address: int, count: int) → List[int]

Read specified number of register starting with given address and return the values from each register in a list.

Parameters

- **address** – address of the first register
- **count** – count of registers to read

Returns list of `int` values

write_registers(*address*: int, *values*: Union[List[int], int])

Write values from the specified address forward.

Parameters

- **address** – address of the first register
- **values** – list with all values

class ModbusTcpCommunicationConfig(*host*: str, *unit*: int, *port*: int = 502)

Bases: object

Configuration dataclass for *ModbusTcpCommunication*.

clean_values()**force_value**(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str

Host is the IP address of the connected device.

is_configdataclass = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 502

TCP port

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

unit: int


Unit number to be used when connecting with Modbus/TCP. Typically this is used when connecting to a relay having Modbus/RTU-connected devices.

exception ModbusTcpConnectionFailedException(*string*=")

Bases: pymodbus.exceptions.ConnectionException

Exception raised when the connection failed.

hvl_ccb.comm.opc

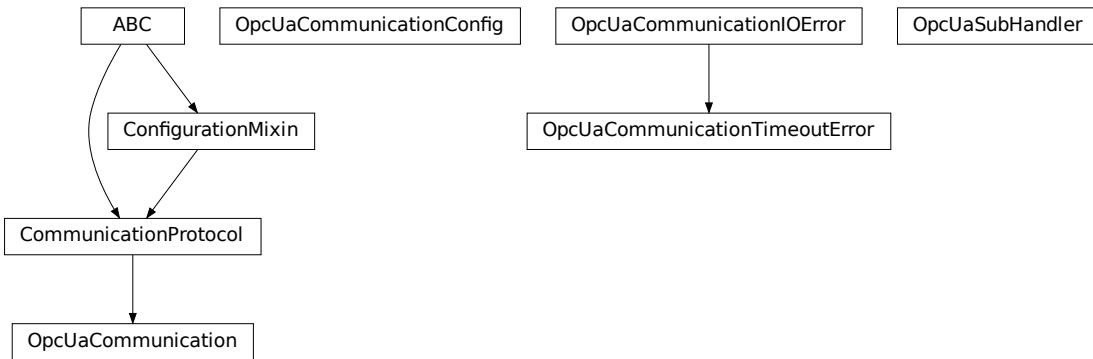

Communication protocol implementing an OPC UA connection. This protocol is used to interface with the “Super-cube” PLC from Siemens.

class OpcUaCommunication(config)

Bases: *hvl_ccb.comm.base.CommunicationProtocol*

Communication protocol implementing an OPC UA connection. Makes use of the package python-opcua.

close() → None

Close the connection to the OPC UA server.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

init_monitored_nodes(node_id: Union[object, Iterable], ns_index: int) → None

Initialize monitored nodes.

Parameters

- **node_id** – one or more strings of node IDs; node IDs are always casted via `str()` method here, hence do not have to be strictly string objects.
- **ns_index** – the namespace index the nodes belong to.

Raises `OpcUaCommunicationIOError` – when protocol was not opened or can't communicate with a OPC UA server

property is_open: bool

Flag indicating if the communication port is open.

Returns `True` if the port is open, otherwise `False`

open() → None

Open the communication to the OPC UA server.

Raises `OpcUaCommunicationIOError` – when communication port cannot be opened.

read(node_id, ns_index)

Read a value from a node with id and namespace index.

Parameters

- **node_id** – the ID of the node to read the value from
- **ns_index** – the namespace index of the node

Returns the value of the node object.

Raises *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

write(*node_id*, *ns_index*, *value*) → None

Write a value to a node with name *name*.

Parameters

- **node_id** – the id of the node to write the value to.
- **ns_index** – the namespace index of the node.
- **value** – the value to write.

Raises *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

```
class OpcUaCommunicationConfig(host: str, endpoint_name: str, port: int = 4840, sub_handler:
                                hvl_ccb.comm.opc.OpcUaSubHandler =
                                <hvl_ccb.comm.opc.OpcUaSubHandler object>, update_period: int = 500,
                                wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int =
                                5)
```

Bases: *object*

Configuration dataclass for OPC UA Communciation.

clean_values()

endpoint_name: str

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str

Hostname or IP-Address of the OPC UA server.

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

max_timeout_retry_nr: int = 5

Maximal number of call re-tries on underlying OPC UA client timeout error

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 4840

Port of the OPC UA server to connect to.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.comm.opc.OpcUaSubHandler object>

object to use for handling subscriptions.

update_period: int = 500

Update period for generating datachange events in OPC UA [milli seconds]

wait_timeout_retry_sec: Union[int, float] = 1

Wait time between re-trying calls on underlying OPC UA client timeout error

exception OpcUaCommunicationIOError

Bases: OSError

OPC-UA communication I/O error.

exception OpcUaCommunicationTimeoutError

Bases: *hvl_ccb.comm.opc.OpcUaCommunicationIOError*

OPC-UA communication timeout error.

class OpcUaSubHandler


Bases: object

Base class for subscription handling of OPC events and data change events. Override methods from this class to add own handling capabilities.

To receive events from server for a subscription data_change and event methods are called directly from receiving thread. Do not do expensive, slow or network operation there. Create another thread if you need to do such a thing.

datachange_notification(node, val, data)

event_notification(event)

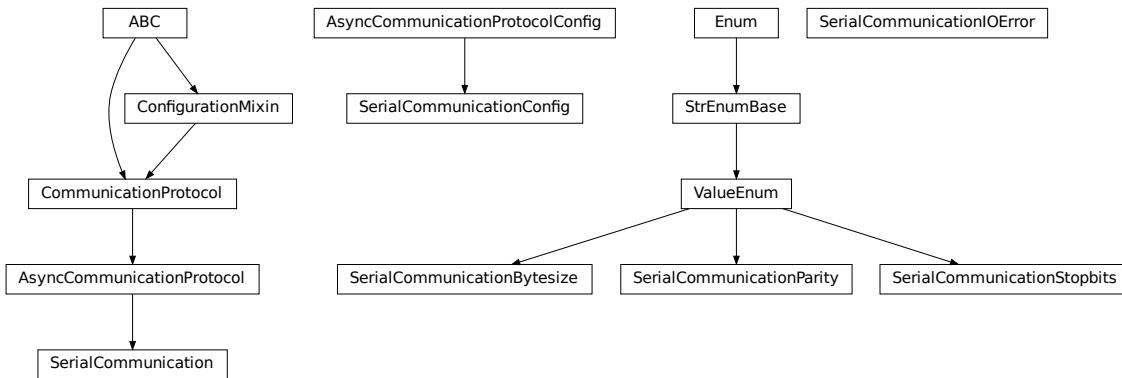hvl_ccb.comm.serial

Communication protocol for serial ports. Makes use of the `pySerial` library.

class SerialCommunication(configuration)
Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocol`

Implements the Communication Protocol for serial ports.

close()

Close the serial connection.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

property is_open: bool

Flag indicating if the serial port is open.

Returns `True` if the serial port is open, otherwise `False`

open()

Open the serial connection.

Raises `SerialCommunicationIOError` – when communication port cannot be opened.

read_bytes() → bytes

Read the bytes from the serial port till the terminator is found. The input buffer may hold additional lines afterwards.

This method uses `self.access_lock` to ensure thread-safety.

Returns Bytes read from the serial port; `b''` if there was nothing to read.

Raises `SerialCommunicationIOError` – when communication port is not opened

read_single_bytes(size: int = 1) → bytes

Read the specified number of bytes from the serial port. The input buffer may hold additional data afterwards.

Returns Bytes read from the serial port; `b''` if there was nothing to read.

write_bytes(data: bytes) → int

Write bytes to the serial port.

This method uses `self.access_lock` to ensure thread-safety.

Parameters `data` – data to write to the serial port

Returns number of bytes written

Raises `SerialCommunicationIOError` – when communication port is not opened

```
class SerialCommunicationBytesize(value=<no_arg>, names=None, module=None, type=None, start=1,  
                                   boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

Serial communication bytesize.

EIGHTBITS = 8

FIVEBITS = 5

SEVENBITS = 7

SIXBITS = 6

```
class SerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',  
                               encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:  
                               Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,  
                               port: Optional[str] = None, baudrate: int = 9600, parity: Union[str,  
                               hvl_ccb.comm.serial.SerialCommunicationParity] =  
                               SerialCommunicationParity.NONE, stopbits: Union[int, float,  
                               hvl_ccb.comm.serial.SerialCommunicationStopbits] =  
                               SerialCommunicationStopbits.ONE, bytesize: Union[int,  
                               hvl_ccb.comm.serial.SerialCommunicationBytesize] =  
                               SerialCommunicationBytesize.EIGHTBITS, timeout: Union[int, float] = 2)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

Configuration dataclass for `SerialCommunication`.

Bytesize

alias of `hvl_ccb.comm.serial.SerialCommunicationBytesize`

Parity

alias of `hvl_ccb.comm.serial.SerialCommunicationParity`

Stopbits

alias of `hvl_ccb.comm.serial.SerialCommunicationStopbits`

baudrate: int = 9600

Baudrate of the serial port

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

Size of a byte, 5 to 8

clean_values()

create_serial_port() → serial.serialposix.Serial

Create a serial port instance according to specification in this configuration

Returns Closed serial port instance

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, *hvl_ccb.comm.serial.SerialCommunicationParity*] = 'N'

Parity to be used for the connection.

port: Optional[str] = None

Port is a string referring to a COM-port (e.g. 'COM3') or a URL. The full list of capabilities is found [on the pyserial documentation](#).

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, float, *hvl_ccb.comm.serial.SerialCommunicationStopbits*] = 1

Stopbits setting, can be 1, 1.5 or 2.

terminator_str() → str

timeout: Union[int, float] = 2

Timeout in seconds for the serial port

exception SerialCommunicationIOError

Bases: OSError

Serial communication related I/O errors.

class SerialCommunicationParity(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

Serial communication parity.

EVEN = 'E'

MARK = 'M'

NAMES = {'E': 'Even', 'M': 'Mark', 'N': 'None', 'O': 'Odd', 'S': 'Space'}

NONE = 'N'

ODD = 'O'

SPACE = 'S'

class SerialCommunicationStopbits(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)


Bases: *hvl_ccb.utils.enum.ValueEnum*

Serial communication stopbits.

ONE = 1

```
ONE_POINT_FIVE = 1.5
TWO = 2
```

[hvl_ccb.comm.tcp](#)

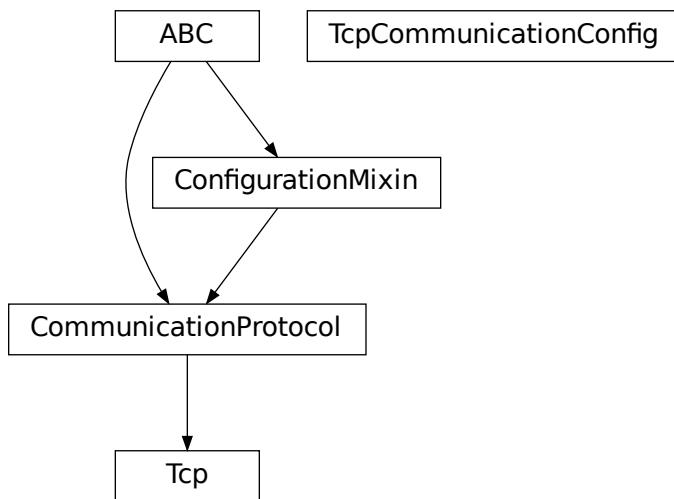

TCP communication protocol.

Makes use of the socket library.

```
class Tcp(configuration)
    Bases: hvl_ccb.comm.base.CommunicationProtocol

    Tcp Communication Protocol.

    close() → None
        Close TCP connection.

    static config_cls() → Type[hvl_ccb.comm.tcp.TcpCommunicationConfig]
        Return the default configdataclass class.

        Returns a reference to the default configdataclass class

    open() → None
        Open TCP connection.

    read() → str
        TCP read function :return: information read from TCP buffer formatted as string

    write(command: str = "") → None
        TCP write function :param command: command string to be sent :return: none

class TcpCommunicationConfig(host: str, port: int = 54321, bufsize: int = 1024)
    Bases: object

    Configuration dataclass for TcpCommunication.

    bufsize: int = 1024
```

clean_values() → None

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.


port: int = 54321

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

hvl_ccb.comm.telnet

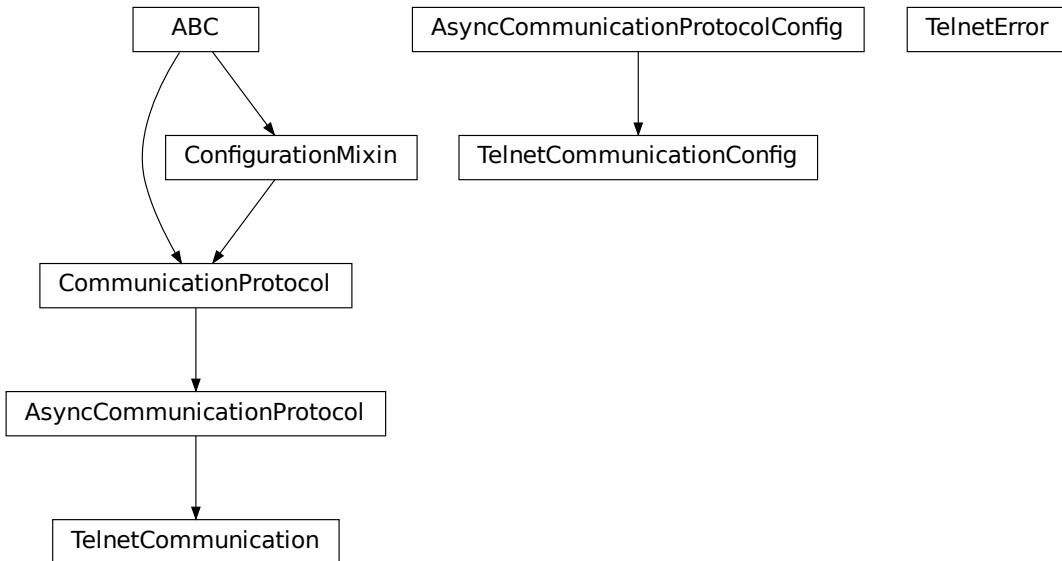

Communication protocol for telnet. Makes use of the `telnetlib` library.

```
class TelnetCommunication(configuration)
    Bases: hvl_ccb.comm.base.AsyncCommunicationProtocol

    Implements the Communication Protocol for telnet.

    close()
        Close the telnet connection unless it is not closed.

    static config_cls()
        Return the default configdataclass class

        Returns a reference to the default configdataclass class

    property is_open: bool
        Is the connection open?

        Returns True for an open connection

    open()
        Open the telnet connection unless it is not yet opened.

    read_bytes() → bytes
        Read data as bytes from the telnet connection.

        Returns data from telnet connection

        Raises TelnetError – when connection is not open, raises an Error during the communication

    write_bytes(data: bytes)
        Write the data as bytes to the telnet connection.

        Parameters data – Data to be sent.
```

Raises `TelnetError` – when connection is not open, raises an Error during the communication

```
class TelnetCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,
                                host: Optional[str] = None, port: int = 0, timeout: Union[int, float] = 0.2)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

Configuration dataclass for `TelnetCommunication`.

clean_values()

create_telnet() → `Optional[telnetlib.Telnet]`

Create a telnet client :return: Opened Telnet object or None if connection is not possible

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: Optional[str] = None

Host to connect to can be localhost or

classmethod keys() → `Sequence[str]`

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → `Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 0

Port at which the host is listening

classmethod required_keys() → `Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

timeout: Union[int, float] = 0.2


Timeout for reading a line

exception TelnetError

Bases: `Exception`

Telnet communication related errors.

hvl_ccb.comm.visa

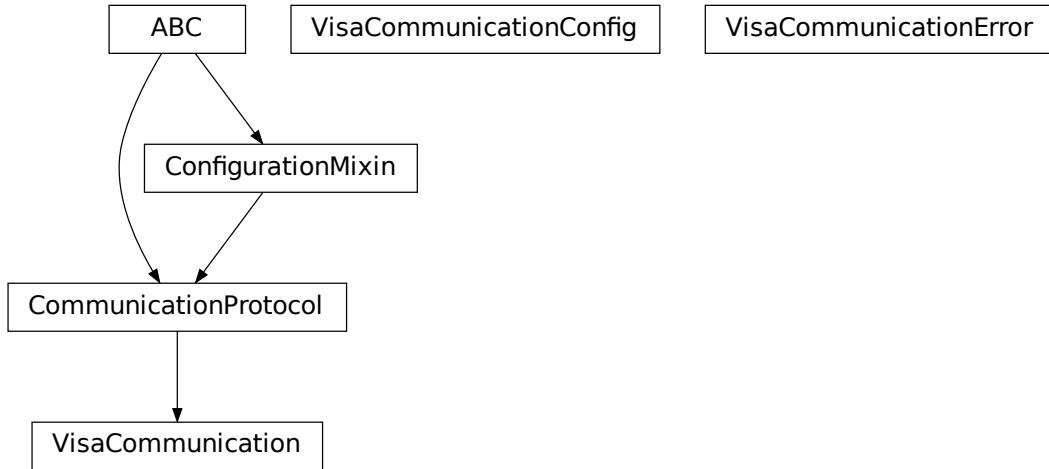

Communication protocol for VISA. Makes use of the pyvisa library. The backend can be NI-Visa or pyvisa-py.

Information on how to install a VISA backend can be found here: https://pyvisa.readthedocs.io/en/master/getting_nivisa.html

So far only TCPIP SOCKET and TCPIP INSTR interfaces are supported.

class VisaCommunication(configuration)

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Implements the Communication Protocol for VISA / SCPI.

MULTI_COMMANDS_MAX = 5

The maximum of commands that can be sent in one round is 5 according to the VISA standard.

MULTI_COMMANDS_SEPARATOR = ';'

The character to separate two commands is ; according to the VISA standard.

WAIT_AFTER_WRITE = 0.08

Small pause in seconds to wait after write operations, allowing devices to really do what we tell them before continuing with further tasks.

close() → None

Close the VISA connection and invalidates the handle.

static config_cls() → Type[`hvl_ccb.comm.visa.VisaCommunicationConfig`]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

open() → None

Open the VISA connection and create the resource.

query(*commands: str) → Union[str, Tuple[str, ...]]

A combination of write(message) and read.

Parameters `commands` – list of commands

Returns list of values

Raises `VisaCommunicationError` – when connection was not started, or when trying to issue too many commands at once.

spoll() → int
 Execute serial poll on the device. Reads the status byte register STB. This is a fast function that can be executed periodically in a polling fashion.

Returns integer representation of the status byte

Raises `VisaCommunicationError` – when connection was not started

write(*commands: str) → None
 Write commands. No answer is read or expected.

Parameters `commands` – one or more commands to send

Raises `VisaCommunicationError` – when connection was not started

```
class VisaCommunicationConfig(host: str, interface_type: Union[str,
    hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType], board: int =
    0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800,
    open_timeout: int = 1000, write_termination: str = '\n', read_termination: str =
    '\n', visa_backend: str = '')
```

Bases: `object`

VisaCommunication configuration dataclass.

```
class InterfaceType(value=<no_arg>, names=None, module=None, type=None, start=1,
    boundary=None)
```

Bases: `hvl_ccb.utils.enum.AutoNumberNameEnum`

Supported VISA Interface types.

TCPIP_INSTR = 2
 VXI-11 protocol

TCPIP_SOCKET = 1
 VISA-Raw protocol

address(host: str, port: Optional[int] = None, board: Optional[int] = None) → str
 Address string specific to the VISA interface type.

Parameters

- `host` – host IP address
- `port` – optional TCP port
- `board` – optional board number

Returns address string

property address: str
 Address string depending on the VISA protocol's configuration.

Returns address string corresponding to current configuration

board: int = 0
 Board number is typically 0 and comes from old bus systems.

chunk_size: int = 204800
 Chunk size is the allocated memory for read operations. The standard is 20kB, and is increased per default here to 200kB. It is specified in bytes.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: str

IP address of the VISA device. DNS names are currently unsupported.

interface_type: Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]

Interface type of the VISA connection, being one of *InterfaceType*.

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

open_timeout: int = 1000

Timeout for opening the connection, in milli seconds.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 5025

TCP port, standard is 5025.

read_termination: str = '\n'

Read termination character.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

timeout: int = 5000

Timeout for commands in milli seconds.

visa_backend: str = ''

Specifies the path to the library to be used with PyVISA as a backend. Defaults to None, which is NI-VISA (if installed), or pyvisa-py (if NI-VISA is not found). To force the use of pyvisa-py, specify '@py' here.

write_termination: str = '\n'

Write termination character.

exception VisaCommunicationError

Bases: Exception

Base class for VisaCommunication errors.

Module contents

Communication protocols subpackage.


`hvl_ccb.dev`

Subpackages

`hvl_ccb.dev.supercube`

Submodules

`hvl_ccb.dev.supercube.base`

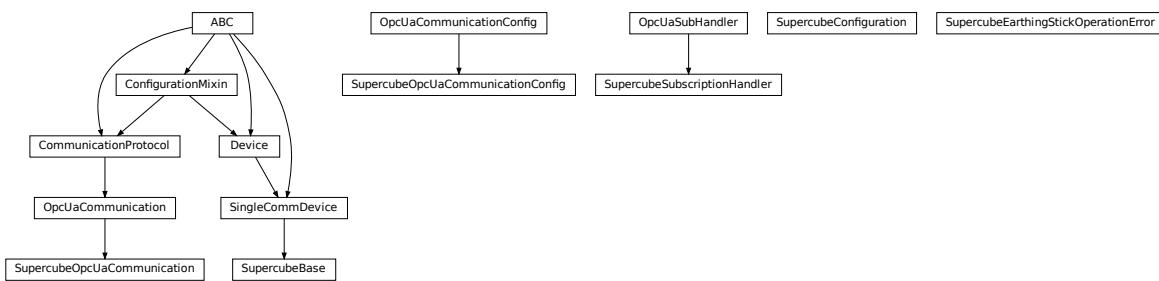

Base classes for the Supercube device.

```
class SupercubeBase(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice
```

Base class for Supercube variants.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

display_message_board() → None

Display 15 newest messages

display_status_board() → None

Display status board.

get_cee16_socket() → bool

Read the on-state of the IEC CEE16 three-phase power socket.

Returns the on-state of the CEE16 power socket

get_door_status(door: int) → `hvl_ccb.dev.supercube.constants.DoorStatus`

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

Parameters `door` – the door number (1..3)

Returns the door status

get_earthing_rod_status(*earthing_rod: int*) → *hvl_ccb.dev.supercube.constants.EarthingRodStatus*

Get the status of a earthing rod. See `constants.EarthingRodStatus` for possible returned earthing rod statuses.

Parameters **earthing_rod** – the earthing rod number (1..3)

Returns the earthing rod status

get_earthing_stick_manual(*number: int*) → *hvl_ccb.dev.supercube.constants.EarthingStickOperation*

Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

Parameters **number** – number of the earthing stick (1..6)

Returns operation of the earthing stick in a manual operating mode (open == 0, close == 1)

Raises `ValueError` – when earthing stick number is not valid

get_earthing_stick_operating_status(*number: int*) →

hvl_ccb.dev.supercube.constants.EarthingStickOperatingStatus

Get the operating status of an earthing stick.

Parameters **number** – number of the earthing stick (1..6)

Returns earthing stick operating status (auto == 0, manual == 1)

Raises `ValueError` – when earthing stick number is not valid

get_earthing_stick_status(*number: int*) → *hvl_ccb.dev.supercube.constants.EarthingStickStatus*

Get the status of an earthing stick, whether it is closed, open or undefined (moving).

Parameters **number** – number of the earthing stick (1..6)

Returns earthing stick status

Raises `ValueError` – when earthing stick number is not valid

get_measurement_ratio(*channel: int*) → float

Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

Parameters **channel** – number of the input channel (1..4)

Returns the ratio

Raises `ValueError` – when channel is not valid

get_measurement_voltage(*channel: int*) → float

Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

Parameters **channel** – number of the input channel (1..4)

Returns measured voltage

Raises `ValueError` – when channel is not valid

get_status() → *hvl_ccb.dev.supercube.constants.SafetyStatus*

Get the safety circuit status of the Supercube. :return: the safety status of the supercube's state machine.

get_support_input(*port: int, contact: int*) → bool

Get the state of a support socket input.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital input read state

Raises ValueError – when port or contact number is not valid

get_support_output(*port: int, contact: int*) → bool

Get the state of a support socket output.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital output read state

Raises ValueError – when port or contact number is not valid

get_t13_socket(*port: int*) → bool

Read the state of a SEV T13 power socket.

Parameters port – is the socket number, one of *constants.T13_SOCKET_PORTS*

Returns on-state of the power socket

Raises ValueError – when port is not valid

operate(*state: bool*) → None

Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

Parameters state – set operate state

operate_earthing_stick(*number: int, operation:*

hvl_ccb.dev.supercube.constants.EarthingStickOperation) → None

Operation of an earthing stick, which is set to manual operation. If an earthing stick is set to manual, it stays closed even if the system is in states RedReady or RedOperate.

Parameters

- **number** – number of the earthing stick (1..6)
- **operation** – earthing stick manual status (close or open)

Raises SupercubeEarthingStickOperationError – when operating status of given number's earthing stick is not manual

quit_error() → None

Quits errors that are active on the Supercube.

read(*node_id: str*)

Local wrapper for the OPC UA communication protocol read method.

Parameters node_id – the id of the node to read.

Returns the value of the variable

ready(*state: bool*) → None

Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

Parameters state – set ready state

set_cee16_socket(*state: bool*) → None

Switch the IEC CEE16 three-phase power socket on or off.

Parameters state – desired on-state of the power socket

Raises ValueError – if state is not of type bool

set_message_board(*msgs: List[str]*, *display_board: bool = True*) → None

Fills messages into message board that display that 15 newest messages with a timestamp.

Parameters

- **msgs** – list of strings
- **display_board** – display 15 newest messages if *True* (default)

Raises ValueError – if there are too many messages or the positions indices are invalid.

set_remote_control(*state: bool*) → None

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

Parameters state – desired remote control state

set_status_board(*msgs: List[str]*, *pos: Optional[List[int]] = None*, *clear_board: bool = True*, *display_board: bool = True*) → None

Sets and displays a status board. The messages and the position of the message can be defined.

Parameters

- **msgs** – list of strings
- **pos** – list of integers [0...14]
- **clear_board** – clear unspecified lines if *True* (default), keep otherwise
- **display_board** – display new status board if *True* (default)

Raises ValueError – if there are too many messages or the positions indices are invalid.

set_support_output(*port: int*, *contact: int*, *state: bool*) → None

Set the state of a support output socket.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **state** – is the desired state of the support output

Raises ValueError – when port or contact number is not valid

set_support_output_impulse(*port: int*, *contact: int*, *duration: float = 0.2*, *pos_pulse: bool = True*) → None

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the *pos_pulse* argument.

This function is blocking.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **duration** – is the length of the impulse in seconds
- **pos_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

Raises ValueError – when port or contact number is not valid

set_t13_socket(*port: int, state: bool*) → None

Set the state of a SEV T13 power socket.

Parameters

- **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*
- **state** – is the desired on-state of the socket

Raises ValueError – when port is not valid or state is not of type bool

start() → None

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

stop() → None

Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

write(*node_id, value*) → None

Local wrapper for the OPC UA communication protocol write method.

Parameters

- **node_id** – the id of the node to read
- **value** – the value to write to the variable

class SupercubeConfiguration(*namespace_index: int = 3, polling_delay_sec: Union[int, float] = 5.0, polling_interval_sec: Union[int, float] = 1.0*)

Bases: object

Configuration dataclass for the Supercube devices.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

namespace_index: int = 3

Namespace of the OPC variables, typically this is 3 (coming from Siemens)

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

polling_delay_sec: Union[int, float] = 5.0

polling_interval_sec: Union[int, float] = 1.0

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception SupercubeEarthingStickOperationError

Bases: Exception

class SupercubeOpcUaCommunication(config)

Bases: [hvl_ccb.comm.opc.OpcUaCommunication](#)

Communication protocol specification for Supercube devices.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class SupercubeOpcUaCommunicationConfig(host: str, endpoint_name: str, port: int = 4840, sub_handler:

*hvl_ccb.comm.opc.OpcUaSubHandler =
<hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler
object>, update_period: int = 500, wait_timeout_retry_sec:
Union[int, float] = 1, max_timeout_retry_nr: int = 5*

Bases: [hvl_ccb.comm.opc.OpcUaCommunicationConfig](#)

Communication protocol configuration for OPC UA, specifications for the Supercube devices.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

**sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
<hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object>**
Subscription handler for data change events

class SupercubeSubscriptionHandler

Bases: [hvl_ccb.comm.opc.OpcUaSubHandler](#)

OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.


datachange_notification(node: opcua.common.node.Node, val, data)

In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.

Parameters

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

hvl_ccb.dev.supercube.constants

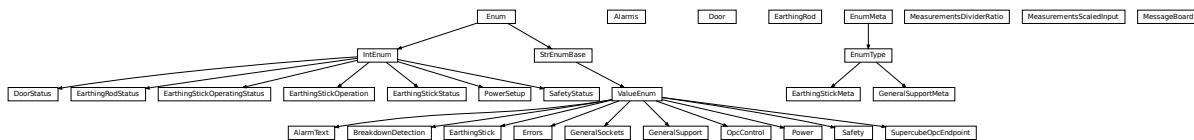

Constants, variable names for the Supercube OPC-connected devices.

```
class AlarmText(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the [AlarmText.get\(\)](#) method to retrieve the enum of an alarm number.

```

Alarm1 = 'STOP Emergency Stop 1'
Alarm10 = 'STOP Earthing stick 2 error while opening'
Alarm11 = 'STOP Earthing stick 3 error while opening'
Alarm12 = 'STOP Earthing stick 4 error while opening'
Alarm13 = 'STOP Earthing stick 5 error while opening'
Alarm14 = 'STOP Earthing stick 6 error while opening'
Alarm15 = 'STOP Earthing stick 1 error while closing'
Alarm16 = 'STOP Earthing stick 2 error while closing'
Alarm17 = 'STOP Earthing stick 3 error while closing'
Alarm18 = 'STOP Earthing stick 4 error while closing'
Alarm19 = 'STOP Earthing stick 5 error while closing'
Alarm2 = 'STOP Emergency Stop 2'
Alarm20 = 'STOP Earthing stick 6 error while closing'
Alarm21 = 'STOP Safety fence 1'
Alarm22 = 'STOP Safety fence 2'
Alarm23 = 'STOP OPC connection error'
Alarm24 = 'STOP Grid power failure'
```

```
Alarm25 = 'STOP UPS failure'
Alarm26 = 'STOP 24V PSU failure'
Alarm3 = 'STOP Emergency Stop 3'
Alarm4 = 'STOP Safety Switch 1 error'
Alarm41 = 'WARNING Door 1: Use earthing rod!'
Alarm42 = 'MESSAGE Door 1: Earthing rod is still in setup.'
Alarm43 = 'WARNING Door 2: Use earthing rod!'
Alarm44 = 'MESSAGE Door 2: Earthing rod is still in setup.'
Alarm45 = 'WARNING Door 3: Use earthing rod!'
Alarm46 = 'MESSAGE Door 3: Earthing rod is still in setup.'
Alarm47 = 'MESSAGE UPS charge < 85%'
Alarm48 = 'MESSAGE UPS running on battery'
Alarm5 = 'STOP Safety Switch 2 error'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error while opening'
```

classmethod get(alarm: int)

Get the attribute of this enum for an alarm number.

Parameters `alarm` – the alarm number

Returns the enum for the desired alarm number

not_defined = 'NO ALARM TEXT DEFINED'

```
class Alarms(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._AlarmEnumBase
```

Alarms enumeration containing all variable NodeID strings for the alarm array.

```
Alarm1 = '"DB_Alarm_HMI"."Alarm1"'
Alarm10 = '"DB_Alarm_HMI"."Alarm10"'
Alarm100 = '"DB_Alarm_HMI"."Alarm100"'
Alarm101 = '"DB_Alarm_HMI"."Alarm101"'
Alarm102 = '"DB_Alarm_HMI"."Alarm102"'
Alarm103 = '"DB_Alarm_HMI"."Alarm103"'
Alarm104 = '"DB_Alarm_HMI"."Alarm104"'
Alarm105 = '"DB_Alarm_HMI"."Alarm105"'
Alarm106 = '"DB_Alarm_HMI"."Alarm106"'
Alarm107 = '"DB_Alarm_HMI"."Alarm107"'
Alarm108 = '"DB_Alarm_HMI"."Alarm108"'
Alarm109 = '"DB_Alarm_HMI"."Alarm109"'
```

```
Alarm11 = '"DB_Alarm_HMI"."Alarm11"'
Alarm110 = '"DB_Alarm_HMI"."Alarm110"'
Alarm111 = '"DB_Alarm_HMI"."Alarm111"'
Alarm112 = '"DB_Alarm_HMI"."Alarm112"'
Alarm113 = '"DB_Alarm_HMI"."Alarm113"'
Alarm114 = '"DB_Alarm_HMI"."Alarm114"'
Alarm115 = '"DB_Alarm_HMI"."Alarm115"'
Alarm116 = '"DB_Alarm_HMI"."Alarm116"'
Alarm117 = '"DB_Alarm_HMI"."Alarm117"'
Alarm118 = '"DB_Alarm_HMI"."Alarm118"'
Alarm119 = '"DB_Alarm_HMI"."Alarm119"'
Alarm12 = '"DB_Alarm_HMI"."Alarm12"'
Alarm120 = '"DB_Alarm_HMI"."Alarm120"'
Alarm121 = '"DB_Alarm_HMI"."Alarm121"'
Alarm122 = '"DB_Alarm_HMI"."Alarm122"'
Alarm123 = '"DB_Alarm_HMI"."Alarm123"'
Alarm124 = '"DB_Alarm_HMI"."Alarm124"'
Alarm125 = '"DB_Alarm_HMI"."Alarm125"'
Alarm126 = '"DB_Alarm_HMI"."Alarm126"'
Alarm127 = '"DB_Alarm_HMI"."Alarm127"'
Alarm128 = '"DB_Alarm_HMI"."Alarm128"'
Alarm129 = '"DB_Alarm_HMI"."Alarm129"'
Alarm13 = '"DB_Alarm_HMI"."Alarm13"'
Alarm130 = '"DB_Alarm_HMI"."Alarm130"'
Alarm131 = '"DB_Alarm_HMI"."Alarm131"'
Alarm132 = '"DB_Alarm_HMI"."Alarm132"'
Alarm133 = '"DB_Alarm_HMI"."Alarm133"'
Alarm134 = '"DB_Alarm_HMI"."Alarm134"'
Alarm135 = '"DB_Alarm_HMI"."Alarm135"'
Alarm136 = '"DB_Alarm_HMI"."Alarm136"'
Alarm137 = '"DB_Alarm_HMI"."Alarm137"'
Alarm138 = '"DB_Alarm_HMI"."Alarm138"'
Alarm139 = '"DB_Alarm_HMI"."Alarm139"'
Alarm14 = '"DB_Alarm_HMI"."Alarm14"'
Alarm140 = '"DB_Alarm_HMI"."Alarm140"'
Alarm141 = '"DB_Alarm_HMI"."Alarm141"'
```

```
Alarm142 = '"DB_Alarm_HMI"."Alarm142"'
Alarm143 = '"DB_Alarm_HMI"."Alarm143"'
Alarm144 = '"DB_Alarm_HMI"."Alarm144"'
Alarm145 = '"DB_Alarm_HMI"."Alarm145"'
Alarm146 = '"DB_Alarm_HMI"."Alarm146"'
Alarm147 = '"DB_Alarm_HMI"."Alarm147"'
Alarm148 = '"DB_Alarm_HMI"."Alarm148"'
Alarm149 = '"DB_Alarm_HMI"."Alarm149"'
Alarm15 = '"DB_Alarm_HMI"."Alarm15"'
Alarm150 = '"DB_Alarm_HMI"."Alarm150"'
Alarm151 = '"DB_Alarm_HMI"."Alarm151"'
Alarm16 = '"DB_Alarm_HMI"."Alarm16"'
Alarm17 = '"DB_Alarm_HMI"."Alarm17"'
Alarm18 = '"DB_Alarm_HMI"."Alarm18"'
Alarm19 = '"DB_Alarm_HMI"."Alarm19"'
Alarm2 = '"DB_Alarm_HMI"."Alarm2"'
Alarm20 = '"DB_Alarm_HMI"."Alarm20"'
Alarm21 = '"DB_Alarm_HMI"."Alarm21"'
Alarm22 = '"DB_Alarm_HMI"."Alarm22"'
Alarm23 = '"DB_Alarm_HMI"."Alarm23"'
Alarm24 = '"DB_Alarm_HMI"."Alarm24"'
Alarm25 = '"DB_Alarm_HMI"."Alarm25"'
Alarm26 = '"DB_Alarm_HMI"."Alarm26"'
Alarm27 = '"DB_Alarm_HMI"."Alarm27"'
Alarm28 = '"DB_Alarm_HMI"."Alarm28"'
Alarm29 = '"DB_Alarm_HMI"."Alarm29"'
Alarm3 = '"DB_Alarm_HMI"."Alarm3"'
Alarm30 = '"DB_Alarm_HMI"."Alarm30"'
Alarm31 = '"DB_Alarm_HMI"."Alarm31"'
Alarm32 = '"DB_Alarm_HMI"."Alarm32"'
Alarm33 = '"DB_Alarm_HMI"."Alarm33"'
Alarm34 = '"DB_Alarm_HMI"."Alarm34"'
Alarm35 = '"DB_Alarm_HMI"."Alarm35"'
Alarm36 = '"DB_Alarm_HMI"."Alarm36"'
Alarm37 = '"DB_Alarm_HMI"."Alarm37"'
Alarm38 = '"DB_Alarm_HMI"."Alarm38"'
```

```
Alarm39 = '"DB_Alarm_HMI"."Alarm39"'
Alarm4 = '"DB_Alarm_HMI"."Alarm4"'
Alarm40 = '"DB_Alarm_HMI"."Alarm40"'
Alarm41 = '"DB_Alarm_HMI"."Alarm41"'
Alarm42 = '"DB_Alarm_HMI"."Alarm42"'
Alarm43 = '"DB_Alarm_HMI"."Alarm43"'
Alarm44 = '"DB_Alarm_HMI"."Alarm44"'
Alarm45 = '"DB_Alarm_HMI"."Alarm45"'
Alarm46 = '"DB_Alarm_HMI"."Alarm46"'
Alarm47 = '"DB_Alarm_HMI"."Alarm47"'
Alarm48 = '"DB_Alarm_HMI"."Alarm48"'
Alarm49 = '"DB_Alarm_HMI"."Alarm49"'
Alarm5 = '"DB_Alarm_HMI"."Alarm5"'
Alarm50 = '"DB_Alarm_HMI"."Alarm50"'
Alarm51 = '"DB_Alarm_HMI"."Alarm51"'
Alarm52 = '"DB_Alarm_HMI"."Alarm52"'
Alarm53 = '"DB_Alarm_HMI"."Alarm53"'
Alarm54 = '"DB_Alarm_HMI"."Alarm54"'
Alarm55 = '"DB_Alarm_HMI"."Alarm55"'
Alarm56 = '"DB_Alarm_HMI"."Alarm56"'
Alarm57 = '"DB_Alarm_HMI"."Alarm57"'
Alarm58 = '"DB_Alarm_HMI"."Alarm58"'
Alarm59 = '"DB_Alarm_HMI"."Alarm59"'
Alarm6 = '"DB_Alarm_HMI"."Alarm6"'
Alarm60 = '"DB_Alarm_HMI"."Alarm60"'
Alarm61 = '"DB_Alarm_HMI"."Alarm61"'
Alarm62 = '"DB_Alarm_HMI"."Alarm62"'
Alarm63 = '"DB_Alarm_HMI"."Alarm63"'
Alarm64 = '"DB_Alarm_HMI"."Alarm64"'
Alarm65 = '"DB_Alarm_HMI"."Alarm65"'
Alarm66 = '"DB_Alarm_HMI"."Alarm66"'
Alarm67 = '"DB_Alarm_HMI"."Alarm67"'
Alarm68 = '"DB_Alarm_HMI"."Alarm68"'
Alarm69 = '"DB_Alarm_HMI"."Alarm69"'
Alarm7 = '"DB_Alarm_HMI"."Alarm7"'
Alarm70 = '"DB_Alarm_HMI"."Alarm70"'
```

```
Alarm71 = '"DB_Alarm_HMI"."Alarm71"'
Alarm72 = '"DB_Alarm_HMI"."Alarm72"'
Alarm73 = '"DB_Alarm_HMI"."Alarm73"'
Alarm74 = '"DB_Alarm_HMI"."Alarm74"'
Alarm75 = '"DB_Alarm_HMI"."Alarm75"'
Alarm76 = '"DB_Alarm_HMI"."Alarm76"'
Alarm77 = '"DB_Alarm_HMI"."Alarm77"'
Alarm78 = '"DB_Alarm_HMI"."Alarm78"'
Alarm79 = '"DB_Alarm_HMI"."Alarm79"'
Alarm8 = '"DB_Alarm_HMI"."Alarm8"'
Alarm80 = '"DB_Alarm_HMI"."Alarm80"'
Alarm81 = '"DB_Alarm_HMI"."Alarm81"'
Alarm82 = '"DB_Alarm_HMI"."Alarm82"'
Alarm83 = '"DB_Alarm_HMI"."Alarm83"'
Alarm84 = '"DB_Alarm_HMI"."Alarm84"'
Alarm85 = '"DB_Alarm_HMI"."Alarm85"'
Alarm86 = '"DB_Alarm_HMI"."Alarm86"'
Alarm87 = '"DB_Alarm_HMI"."Alarm87"'
Alarm88 = '"DB_Alarm_HMI"."Alarm88"'
Alarm89 = '"DB_Alarm_HMI"."Alarm89"'
Alarm9 = '"DB_Alarm_HMI"."Alarm9"'
Alarm90 = '"DB_Alarm_HMI"."Alarm90"'
Alarm91 = '"DB_Alarm_HMI"."Alarm91"'
Alarm92 = '"DB_Alarm_HMI"."Alarm92"'
Alarm93 = '"DB_Alarm_HMI"."Alarm93"'
Alarm94 = '"DB_Alarm_HMI"."Alarm94"'
Alarm95 = '"DB_Alarm_HMI"."Alarm95"'
Alarm96 = '"DB_Alarm_HMI"."Alarm96"'
Alarm97 = '"DB_Alarm_HMI"."Alarm97"'
Alarm98 = '"DB_Alarm_HMI"."Alarm98"'
Alarm99 = '"DB_Alarm_HMI"."Alarm99"'

class BreakdownDetection(value=<no_arg>, names=None, module=None, type=None, start=1,
                        boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Node ID strings for the breakdown detection.

TODO: these variable NodeIDs are not tested and/or correct yet.
```

```
activated = '"Ix_Allg_Breakdown_activated"
Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the system or not.

reset = '"Qx_Allg_Breakdown_reset"
Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

triggered = '"Ix_Allg_Breakdown_triggered"
Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using the safety circuit state, therefore no method is implemented to read this out directly.

class Door(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._DoorEnumBase

Variable NodeID strings for doors.

status_1 = '"DB_Safety_Circuit"."Door_1"."si_HMI_status"
status_2 = '"DB_Safety_Circuit"."Door_2"."si_HMI_status"
status_3 = '"DB_Safety_Circuit"."Door_3"."si_HMI_status"

class DoorStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for doors.

closed = 2
Door is closed, but not locked.

error = 4
Door has an error or was opened in locked state (either with emergency stop or from the inside).

inactive = 0
not enabled in Supercube HMI setup, this door is not supervised.

locked = 3
Door is closed and locked (safe state).

open = 1
Door is open.

class EarthingRod(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._DoorEnumBase

Variable NodeID strings for earthing rods.

status_1 = '"DB_Safety_Circuit"."Door_1"."Ix_earthingrod"
status_2 = '"DB_Safety_Circuit"."Door_2"."Ix_earthingrod"
status_3 = '"DB_Safety_Circuit"."Door_3"."Ix_earthingrod"

class EarthingRodStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: aenum.IntEnum

Possible status values for earthing rods.

experiment_blocked = 0
earthing rod is somewhere in the experiment and blocks the start of the experiment

experiment_ready = 1
earthing rod is hanging next to the door, experiment is ready to operate
```

```
class EarthingStick(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the earthing in manual mode.

```
classmethod manual(number: int)
```

Get the manual enum instance for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the manual instance

Raises **ValueError** – when earthing stick number is not valid

```
manual_1 = '"DB_Safety_Circuit"."Earthstick_1"."sx_earthing_manually"'
```

```
manual_2 = '"DB_Safety_Circuit"."Earthstick_2"."sx_earthing_manually"'
```

```
manual_3 = '"DB_Safety_Circuit"."Earthstick_3"."sx_earthing_manually"'
```

```
manual_4 = '"DB_Safety_Circuit"."Earthstick_4"."sx_earthing_manually"'
```

```
manual_5 = '"DB_Safety_Circuit"."Earthstick_5"."sx_earthing_manually"'
```

```
manual_6 = '"DB_Safety_Circuit"."Earthstick_6"."sx_earthing_manually"'
```

```
classmethod manuals() → Tuple[hvl_ccb.dev.supercube.constants.EarthingStick, ...]
```

Get all earthing stick manual instances.

Returns tuple of manual instances

```
property number: int
```

Get corresponding earthing stick number.

Returns earthing stick number (1..6)

```
classmethod operating_status(number: int)
```

Get the operating status enum instance for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the operating status instance

Raises **ValueError** – when earthing stick number is not valid

```
operating_status_1 = '"DB_Safety_Circuit"."Earthstick_1"."sx_manual_control_active"'
```

```
operating_status_2 = '"DB_Safety_Circuit"."Earthstick_2"."sx_manual_control_active"'
```

```
operating_status_3 = '"DB_Safety_Circuit"."Earthstick_3"."sx_manual_control_active"'
```

```
operating_status_4 = '"DB_Safety_Circuit"."Earthstick_4"."sx_manual_control_active"'
```

```
operating_status_5 = '"DB_Safety_Circuit"."Earthstick_5"."sx_manual_control_active"'
```

```
operating_status_6 = '"DB_Safety_Circuit"."Earthstick_6"."sx_manual_control_active"'
```

```
classmethod operating_statuses() → Tuple[hvl_ccb.dev.supercube.constants.EarthingStick, ...]
```

Get all earthing stick operating status instances.

Returns tuple of operating status instances

```
classmethod range() → Sequence[int]
```

Integer range of all earthing sticks.

Returns sequence of earthing sticks numbers

```
classmethod status(number: int) → hvl_ccb.dev.supercube.constants.EarthingStick
```

Get the status enum instance for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status instance

Raises **ValueError** – when earthing stick number is not valid

```
status_1 = '"DB_Safety_Circuit"."Earthstick_1"."si_HMI_Status"'
```

```
status_2 = '"DB_Safety_Circuit"."Earthstick_2"."si_HMI_Status"'
```

```
status_3 = '"DB_Safety_Circuit"."Earthstick_3"."si_HMI_Status"'
```

```
status_4 = '"DB_Safety_Circuit"."Earthstick_4"."si_HMI_Status"'
```

```
status_5 = '"DB_Safety_Circuit"."Earthstick_5"."si_HMI_Status"'
```

```
status_6 = '"DB_Safety_Circuit"."Earthstick_6"."si_HMI_Status"'
```

```
classmethod statuses() → Tuple[hvl_ccb.dev.supercube.constants.EarthingStick, ...]
```

Get all earthing stick status instances.

Returns tuple of status instances

```
class EarthingStickMeta(clsname, bases, clsdict, **kwargs)
```

Bases: aenum.EnumType

```
class EarthingStickOperatingStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: aenum.IntEnum

Operating Status for an earthing stick. Stick can be used in auto or manual mode.

```
auto = 0
```

```
manual = 1
```

```
class EarthingStickOperation(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: aenum.IntEnum

Operation of the earthing stick in manual operating mode. Can be closed or opened.

```
close = 1
```

```
open = 0
```

```
class EarthingStickStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: aenum.IntEnum

Status of an earthing stick. These are the possible values in the status integer e.g. in *EarthingStick.status_1*.

```
closed = 1
```

Earthing is closed (safe).

```
error = 3
```

Earthing is in error, e.g. when the stick did not close correctly or could not open.

```
inactive = 0
```

Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be enabled in the Supercube HMI setup.

open = 2

Earthing is open (not safe).

class Errors(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

Variable NodeID strings for information regarding error, warning and message handling.

message = "DB_Message_Buffer". "Info_active"

Boolean read-only variable telling if a message is active.

quit = "DB_Message_Buffer". "Reset_button"

Writable boolean for the error quit button.

stop = "DB_Message_Buffer". "Stop_active"

Boolean read-only variable telling if a stop is active.

warning = "DB_Message_Buffer". "Warning_active"

Boolean read-only variable telling if a warning is active.

class GeneralSockets(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

NodeID strings for the power sockets (3x T13 and 1xCEE16).

cee16 = "Qx_Allg_Socket_CEE16"

CEE16 socket (writeable boolean).

t13_1 = "Qx_Allg_Socket_T13_1"

SEV T13 socket No. 1 (writable boolean).

t13_2 = "Qx_Allg_Socket_T13_2"

SEV T13 socket No. 2 (writable boolean).

t13_3 = "Qx_Allg_Socket_T13_3"

SEV T13 socket No. 3 (writable boolean).

class GeneralSupport(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

NodeID strings for the support inputs and outputs.

classmethod contact_range() → Sequence[int]

Integer range of all contacts.

Returns sequence of contact numbers

in_1_1 = "Ix_Allg_Support1_1"

in_1_2 = "Ix_Allg_Support1_2"

in_2_1 = "Ix_Allg_Support2_1"

in_2_2 = "Ix_Allg_Support2_2"

in_3_1 = "Ix_Allg_Support3_1"

in_3_2 = "Ix_Allg_Support3_2"

in_4_1 = "Ix_Allg_Support4_1"

in_4_2 = "Ix_Allg_Support4_2"

in_5_1 = "Ix_Allg_Support5_1"

in_5_2 = "Ix_Allg_Support5_2"

```
in_6_1 = '"Ix_Allg_Support6_1"'
in_6_2 = '"Ix_Allg_Support6_2"'

```

classmethod input(*port: int, contact: int*)
Get the NodeID string for a support input.

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string**Raises ValueError** – when port or contact number is not valid

```
out_1_1 = '"Qx_Allg_Support1_1"'
out_1_2 = '"Qx_Allg_Support1_2"'
out_2_1 = '"Qx_Allg_Support2_1"'
out_2_2 = '"Qx_Allg_Support2_2"'
out_3_1 = '"Qx_Allg_Support3_1"'
out_3_2 = '"Qx_Allg_Support3_2"'
out_4_1 = '"Qx_Allg_Support4_1"'
out_4_2 = '"Qx_Allg_Support4_2"'
out_5_1 = '"Qx_Allg_Support5_1"'
out_5_2 = '"Qx_Allg_Support5_2"'
out_6_1 = '"Qx_Allg_Support6_1"'
out_6_2 = '"Qx_Allg_Support6_2"'

```

classmethod output(*port: int, contact: int*)

Get the NodeID string for a support output.

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string**Raises ValueError** – when port or contact number is not valid**classmethod port_range()** → Sequence[int]

Integer range of all ports.

Returns sequence of port numbers**class GeneralSupportMeta**(*clsname, bases, clsdict, **kwargs*)

Bases: aenum.EnumType

class MeasurementsDividerRatio(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: hvl_ccb.dev.supercube.constants._InputEnumBase

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

```
input_1 = '"DB_Measurements"."si_Divider_Ratio_1"'
input_2 = '"DB_Measurements"."si_Divider_Ratio_2"'
input_3 = '"DB_Measurements"."si_Divider_Ratio_3"'
input_4 = '"DB_Measurements"."si_Divider_Ratio_4"'

class MeasurementsScaledInput(value=<no_arg>, names=None, module=None, type=None, start=1,
                                boundary=None)
Bases: hvl_ccb.dev.supercube.constants._InputEnumBase
```

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in [MeasurementsDividerRatio](#).

```
input_1 = '"DB_Measurements"."si_scaled_Voltage_Input_1"'
input_2 = '"DB_Measurements"."si_scaled_Voltage_Input_2"'
input_3 = '"DB_Measurements"."si_scaled_Voltage_Input_3"'
input_4 = '"DB_Measurements"."si_scaled_Voltage_Input_4"'

class MessageBoard(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._LineEnumBase
```

Variable NodeID strings for message board lines.

```
line_1 = '"DB_OP_C_Connection"."Is_status_Line_1"'
line_10 = '"DB_OP_C_Connection"."Is_status_Line_10"'
line_11 = '"DB_OP_C_Connection"."Is_status_Line_11"'
line_12 = '"DB_OP_C_Connection"."Is_status_Line_12"'
line_13 = '"DB_OP_C_Connection"."Is_status_Line_13"'
line_14 = '"DB_OP_C_Connection"."Is_status_Line_14"'
line_15 = '"DB_OP_C_Connection"."Is_status_Line_15"'
line_2 = '"DB_OP_C_Connection"."Is_status_Line_2"'
line_3 = '"DB_OP_C_Connection"."Is_status_Line_3"'
line_4 = '"DB_OP_C_Connection"."Is_status_Line_4"'
line_5 = '"DB_OP_C_Connection"."Is_status_Line_5"'
line_6 = '"DB_OP_C_Connection"."Is_status_Line_6"'
line_7 = '"DB_OP_C_Connection"."Is_status_Line_7"'
line_8 = '"DB_OP_C_Connection"."Is_status_Line_8"'
line_9 = '"DB_OP_C_Connection"."Is_status_Line_9"'
```

```
class OpcControl(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

Variable NodeID strings for supervision of the OPC connection from the controlling workstation to the Super-cube.

```
active = '"DB_OP_C_Connection"."sx_OP_C_active"'
        writable boolean to enable OPC remote control and display a message window on the Supercube HMI.

live = '"DB_OP_C_Connection"."sx_OP_C_lifebit"'
```

```
class Power(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Variable NodeID strings concerning power data.

TODO: these variable NodeIDs are not tested and/or correct yet, they don't exist yet on Supercube side.

current_primary = 'Qr_Power_FU_actual_Current'
Primary current in ampere, measured by the frequency converter. (read-only)

frequency = 'Ir_Power_FU_Frequency'
Frequency converter output frequency. (read-only)

setup = 'Qi_Power_Setup'
Power setup that is configured using the Supercube HMI. The value corresponds to the ones in PowerSetup. (read-only)

voltage_max = 'Iw_Power_max_Voltage'
Maximum voltage allowed by the current experimental setup. (read-only)

voltage_primary = 'Qr_Power_FU_actual_Voltage'
Primary voltage in volts, measured by the frequency converter at its output. (read-only)

voltage_slope = 'Ir_Power_dUdt'
Voltage slope in V/s.

voltage_target = 'Ir_Power_Target_Voltage'
Target voltage setpoint in V.

class PowerSetup(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible power setups corresponding to the value of variable Power.setup.

AC_DoubleStage_150kV = 4
AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum voltage of 150kV.

AC_DoubleStage_200kV = 5
AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

AC_SingleStage_100kV = 3
AC voltage with MWB transformer set to 100kV maximum voltage.

AC_SingleStage_50kV = 2
AC voltage with MWB transformer set to 50kV maximum voltage.

DC_DoubleStage_280kV = 8
DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single stage transformer with Greinacher voltage doubling rectifier)

DC_SingleStage_140kV = 7
DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

External = 1
External power supply fed through blue CEE32 input using isolation transformer and safety switches of the Supercube, or using an external safety switch attached to the Supercube Type B.

Internal = 6
Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

NoPower = 0
No safety switches, use only safety components (doors, fence, earthing...) without any power.
```

```
class Safety(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

status = '"DB_Safety_Circuit"."si_safe_status"'
Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in SafetyStatus.

switch_to_operate = '"DB_Safety_Circuit"."sx_safe_switch_to_operate"'
Writable boolean for switching to Red Operate (locket, HV on) state.

switch_to_ready = '"DB_Safety_Circuit"."sx_safe_switch_to_ready"'
Writable boolean for switching to Red Ready (locked, HV off) state.

class SafetyStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Safety status values that are possible states returned from hvl_ccb.dev.supercube.base.Supercube.get_status(). These values correspond to the states of the Supercube’s safety circuit statemachine.

Error = 6
System is in error mode.

GreenNotReady = 1
System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

GreenReady = 2
System is safe and all safety elements are in place to be able to switch to ready.

Initializing = 0
System is initializing or booting.

QuickStop = 5
Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

RedOperate = 4
System is locked in red state and in operate mode, i.e. high voltage on.


RedReady = 3
System is locked in red state and ready to go to operate mode.

class SupercubeOpcEndpoint(value=<no_arg>, names=None, module=None, type=None, start=1,
                             boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

OPC Server Endpoint strings for the supercube variants.

A = 'Supercube Typ A'
B = 'Supercube Typ B'

T13_SOCKET_PORTS = (1, 2, 3)
Port numbers of SEV T13 power socket
```

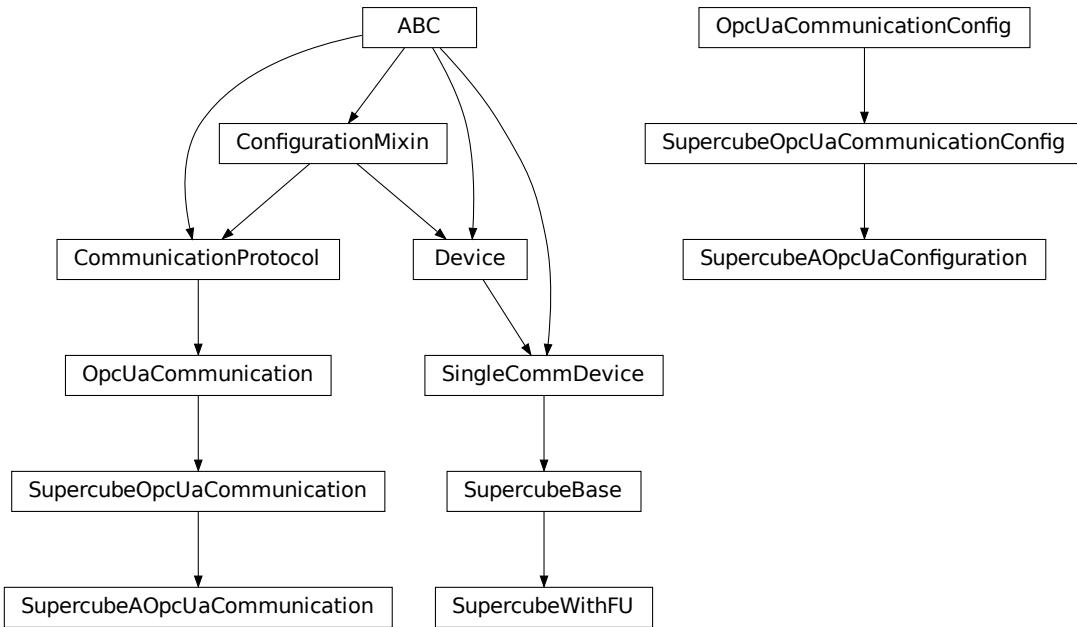hvl_ccb.dev.supercube.typ_a

Supercube Typ A module.

class SupercubeAOpcUaCommunication(config)
Bases: *hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunication*

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class SupercubeAOpcUaConfiguration(host: str, endpoint_name: str = 'Supercube Typ A', port: int = 4840, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object at 0x7fed1b3dd210>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)
Bases: *hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunicationConfig*

endpoint_name: str = 'Supercube Typ A'

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

class SupercubeWithFU(*com, dev_config=None*)
Bases: *hvl_ccb.dev.supercube.base.SupercubeBase*

Variant A of the Supercube with frequency converter.

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

fso_reset() → None
TODO: test fso_reset with device
Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.

get_frequency() → float
TODO: test get_frequency with device
Read the electrical frequency of the current Supercube setup.

Returns the frequency in Hz

get_fso_active() → bool
TODO: test get_fso_active with device
Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.

Returns state of the FSO functionality

get_max_voltage() → float
TODO: test get_max_voltage with device
Reads the maximum voltage of the setup and returns in V.

Returns the maximum voltage of the setup in V.

get_power_setup() → *hvl_ccb.dev.supercube.constants.PowerSetup*
TODO: test get_power_setup with device
Return the power setup selected in the Supercube's settings.

Returns the power setup

get_primary_current() → float
TODO: get_primary_current with device
Read the current primary current at the output of the frequency converter (before transformer).

Returns primary current in A

`get_primary_voltage() → float`

TODO: test `get_primary_voltage` with device

Read the current primary voltage at the output of the frequency converter (before transformer).

Returns primary voltage in V

`get_target_voltage() → float`

TODO: test `get_target_voltage` with device

Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to `set_target_voltage()`.

Returns the setpoint voltage in V.

`set_slope(slope: float) → None`

TODO: test `set_slope` with device

Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.

Parameters `slope` – voltage slope in V/s (0..15)


`set_target_voltage(volt_v: float) → None`

TODO: test `set_target_voltage` with device

Set the output voltage to a defined value in V.

Parameters `volt_v` – the desired voltage in V

hvl_ccb.dev.supercube.typ_b

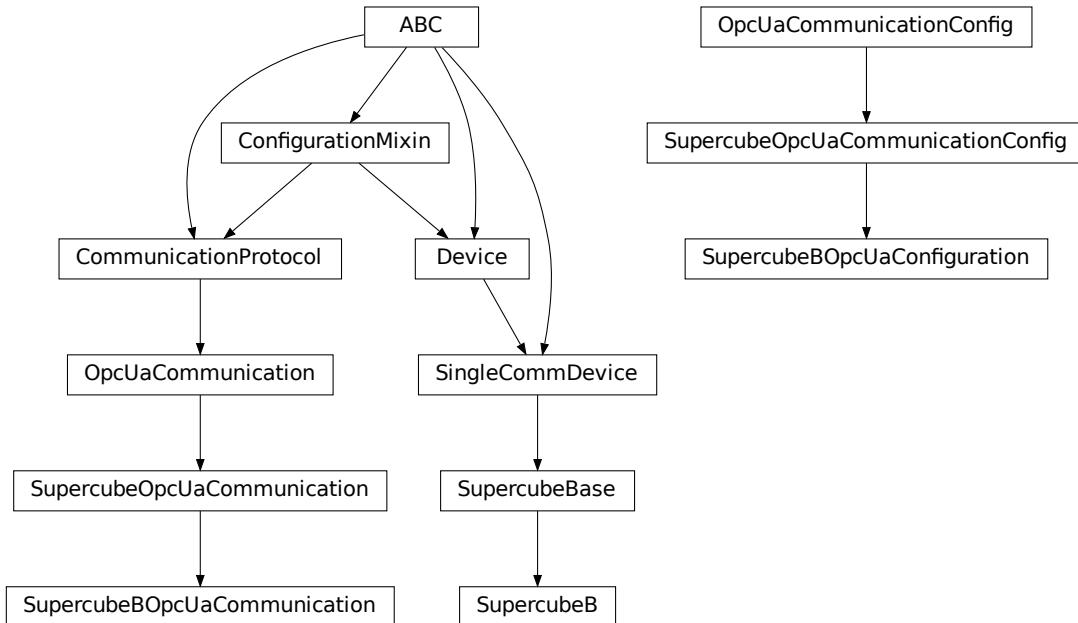

Supercube Typ B module.

```
class SupercubeB(com, dev_config=None)
Bases: hvl_ccb.dev.supercube.base.SupercubeBase

Variant B of the Supercube without frequency converter but external safety switches.

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

class SupercubeB0pcUaCommunication(config)
Bases: hvl_ccb.dev.supercube.base.Supercube0pcUaCommunication

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

class SupercubeB0pcUaConfiguration(host: str, endpoint_name: str = 'Supercube Typ B', port: int = 4840,
                                   sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
                                   <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object at
                                   0x7fed1b3dd210>, update_period: int = 500, wait_timeout_retry_sec:
                                   Union[int, float] = 1, max_timeout_retry_nr: int = 5)
Bases: hvl_ccb.dev.supercube.base.Supercube0pcUaCommunicationConfig

endpoint_name: str = 'Supercube Typ B'
    Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing
    after value has been forced on fieldname.

    Parameters
        • fieldname – name of the field
        • value – value to assign

classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
    on instantiation.

    Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
    instantiation.

    Returns a list of strings containing all required keys.
```


Module contents

Supercube package with implementation for system versions from 2019 on (new concept with hard-PLC Siemens S7-1500 as CPU).

`hvl_ccb.dev.supercube2015`

Submodules

`hvl_ccb.dev.supercube2015.base`

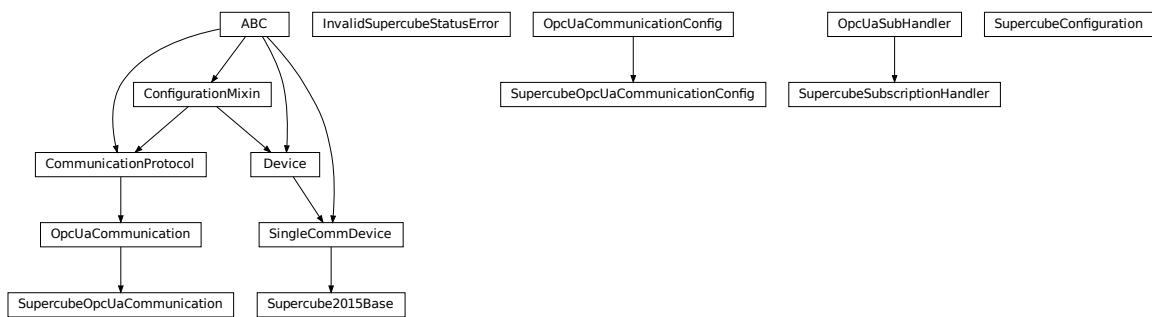

Base classes for the Supercube device.

`exception InvalidSupercubeStatusError`

Bases: `Exception`

Exception raised when supercube has invalid status.

`class Supercube2015Base(com, dev_config=None)`

Bases: `hvl_ccb.dev.base.SingleCommDevice`

Base class for Supercube variants.

`static config_cls()`

Return the default configdataclass class.

Returns a reference to the default configdataclass class

`static default_com_cls()`

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

`get_cee16_socket() → bool`

Read the on-state of the IEC CEE16 three-phase power socket.

Returns the on-state of the CEE16 power socket

`get_door_status(door: int) → hvl_ccb.dev.supercube2015.constants.DoorStatus`

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

Parameters `door` – the door number (1..3)

Returns the door status

get_earthing_manual(*number: int*) → bool

Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

Parameters **number** – number of the earthing stick (1..6)

Returns earthing stick manual status

get_earthing_status(*number: int*) → int

Get the status of an earthing stick, whether it is closed, open or undefined (moving).

Parameters **number** – number of the earthing stick (1..6)

Returns earthing stick status; see constants.EarthlingStickStatus

get_measurement_ratio(*channel: int*) → float

Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

Attention: Supercube 2015 does not have a separate ratio for every analog input. Therefore there is only one ratio for channel = 1.

Parameters **channel** – number of the input channel (1..4)

Returns the ratio

get_measurement_voltage(*channel: int*) → float

Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

Attention: In contrast to the *new* Supercube, the old one returns here the input voltage read at the ADC. It is not scaled by a factor.

Parameters **channel** – number of the input channel (1..4)

Returns measured voltage

get_status() → int

Get the safety circuit status of the Supercube.

Returns the safety status of the supercube's state machine; see *constants.SafetyStatus*.

get_support_input(*port: int, contact: int*) → bool

Get the state of a support socket input.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital input read state

get_support_output(*port: int, contact: int*) → bool

Get the state of a support socket output.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital output read state

get_t13_socket(*port: int*) → bool

Read the state of a SEV T13 power socket.

Parameters **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*

Returns on-state of the power socket

horn(*state: bool*) → None

Turns acoustic horn on or off.

Parameters **state** – Turns horn on (True) or off (False)

operate(*state: bool*) → None

Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

Parameters **state** – set operate state

quit_error() → None

Quits errors that are active on the Supercube.

read(*node_id: str*)

Local wrapper for the OPC UA communication protocol read method.

Parameters **node_id** – the id of the node to read.

Returns the value of the variable

ready(*state: bool*) → None

Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

Parameters **state** – set ready state

set_cee16_socket(*state: bool*) → None

Switch the IEC CEE16 three-phase power socket on or off.

Parameters **state** – desired on-state of the power socket

Raises **ValueError** – if state is not of type bool

set_earthing_manual(*number: int, manual: bool*) → None

Set the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

Parameters

- **number** – number of the earthing stick (1..6)
- **manual** – earthing stick manual status (True or False)

set_remote_control(*state: bool*) → None

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

Parameters **state** – desired remote control state

set_support_output(*port: int, contact: int, state: bool*) → None

Set the state of a support output socket.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **state** – is the desired state of the support output

set_support_output_impulse(*port: int, contact: int, duration: float = 0.2, pos_pulse: bool = True*) → None

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the pos_pulse argument.

This function is blocking.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **duration** – is the length of the impulse in seconds
- **pos_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

set_t13_socket(*port: int, state: bool*) → None

Set the state of a SEV T13 power socket.

Parameters

- **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*
- **state** – is the desired on-state of the socket

start() → None

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

stop() → None

Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

write(*node_id, value*) → None

Local wrapper for the OPC UA communication protocol write method.

Parameters

- **node_id** – the id of the node to read
- **value** – the value to write to the variable

class SupercubeConfiguration(namespace_index: int = 7)

Bases: *object*

Configuration dataclass for the Supercube devices.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

namespace_index: int = 7

Namespace of the OPC variables, typically this is 3 (coming from Siemens)

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

class SupercubeOpcUaCommunication(config)

Bases: *hvl_ccb.comm.opc.OpcUaCommunication*

Communication protocol specification for Supercube devices.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class SupercubeOpcUaCommunicationConfig(host: str, endpoint_name: str, port: int = 4845, sub_handler:

```
hvl_ccb.comm.opc.OpcUaSubHandler =
<hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler
object>, update_period: int = 500, wait_timeout_retry_sec:
Union[int, float] = 1, max_timeout_retry_nr: int = 5
```

Bases: *hvl_ccb.comm.opc.OpcUaCommunicationConfig*

Communication protocol configuration for OPC UA, specifications for the Supercube devices.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 4845

Port of the OPC UA server to connect to.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

```
sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
<hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object>
    Subscription handler for data change events

class SupercubeSubscriptionHandler
Bases: hvl_ccb.comm.opc.OpcUaSubHandler


OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.

datachange_notification(node: opcua.common.node.Node, val, data)
    In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.
```

Parameters

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

hvl_ccb.dev.supercube2015.constants

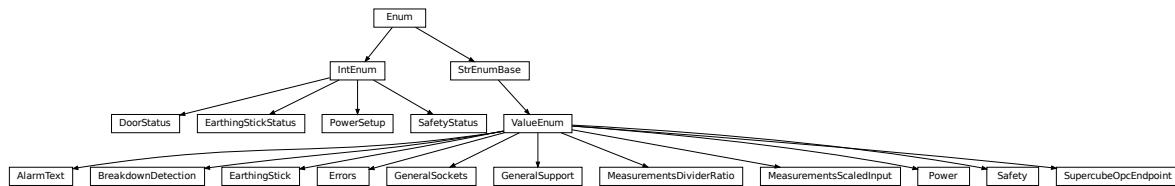

Constants, variable names for the Supercube OPC-connected devices.

```
class AlarmText(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the [AlarmText.get\(\)](#) method to retrieve the enum of an alarm number.

```
Alarm0 = 'No Alarm.'
Alarm1 = 'STOP Safety switch 1 error'
Alarm10 = 'STOP Earthing stick 2 error'
Alarm11 = 'STOP Earthing stick 3 error'
Alarm12 = 'STOP Earthing stick 4 error'
Alarm13 = 'STOP Earthing stick 5 error'
Alarm14 = 'STOP Earthing stick 6 error'
Alarm17 = 'STOP Source switch error'
Alarm19 = 'STOP Fence 1 error'
Alarm2 = 'STOP Safety switch 2 error'
Alarm20 = 'STOP Fence 2 error'
Alarm21 = 'STOP Control error'
```

```

Alarm22 = 'STOP Power outage'
Alarm3 = 'STOP Emergency Stop 1'
Alarm4 = 'STOP Emergency Stop 2'
Alarm5 = 'STOP Emergency Stop 3'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error'

classmethod get(alarm: int)
    Get the attribute of this enum for an alarm number.

        Parameters alarm – the alarm number

        Returns the enum for the desired alarm number

not_defined = 'NO ALARM TEXT DEFINED'

class BreakdownDetection(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Node ID strings for the breakdown detection.

activated = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.Breakdowndetection.connect'
Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the system or not.

reset = 'hvl-ipc.WINAC.Support60utA'
Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

triggered = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.Breakdowndetection.triggered'
Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using the safety circuit state, therefore no method is implemented to read this out directly.

class DoorStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for doors.

closed = 2
Door is closed, but not locked.

error = 4
Door has an error or was opened in locked state (either with emergency stop or from the inside).

inactive = 0
not enabled in Supercube HMI setup, this door is not supervised.

locked = 3
Door is closed and locked (safe state).

open = 1
Door is open.

class EarthingStick(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

```

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the earthing in manual mode.

classmethod manual(*number: int*)

Get the manual enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the manual enum

```
manual_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.MANUAL'  
manual_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.MANUAL'  
manual_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.MANUAL'  
manual_4 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.MANUAL'  
manual_5 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.MANUAL'  
manual_6 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.MANUAL'  
  
status_1_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.CLOSE'  
status_1_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.CONNECT'  
status_1_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.OPEN'  
  
status_2_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.CLOSE'  
status_2_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.CONNECT'  
status_2_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.OPEN'  
  
status_3_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.CLOSE'  
status_3_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.CONNECT'  
status_3_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.OPEN'  
  
status_4_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.CLOSE'  
status_4_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.CONNECT'  
status_4_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.OPEN'  
  
status_5_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.CLOSE'  
status_5_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.CONNECT'  
status_5_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.OPEN'  
  
status_6_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.CLOSE'  
status_6_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.CONNECT'  
status_6_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.OPEN'
```

classmethod status_closed(*number: int*)

Get the status enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status enum

classmethod status_connected(*number: int*)

Get the status enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status enum

classmethod **status_open**(*number*: int)

Get the status enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status enum

class **EarthingStickStatus**(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: aenum.IntEnum

Status of an earthing stick. These are the possible values in the status integer e.g. in EarthingStick.status_1.

closed = 1

Earthing is closed (safe).

error = 3

Earthing is in error, e.g. when the stick did not close correctly or could not open.

inactive = 0

Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be enabled in the Supercube HMI setup.

open = 2

Earthing is open (not safe).

class **Errors**(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: hvl_ccb.utils.enum.ValueEnum

Variable NodeID strings for information regarding error, warning and message handling.

quit = 'hvl-ipc.WINAC.SYSTEMSTATE.Faultconfirmation'

Writable boolean for the error quit button.

stop = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'

Boolean read-only variable telling if a stop is active.

stop_number = 'hvl-ipc.WINAC.SYSTEMSTATE.Errornumber'

class **GeneralSockets**(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: hvl_ccb.utils.enum.ValueEnum

NodeID strings for the power sockets (3x T13 and 1xCEE16).

cee16 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.CEE16'

CEE16 socket (writeable boolean).

t13_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_1'

SEV T13 socket No. 1 (writable boolean).

t13_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_2'

SEV T13 socket No. 2 (writable boolean).

t13_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_3'

SEV T13 socket No. 3 (writable boolean).

class **GeneralSupport**(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: hvl_ccb.utils.enum.ValueEnum

NodeID strings for the support inputs and outputs.

in_1_1 = 'hvl-ipc.WINAC.Support1InA'

```
in_1_2 = 'hvl-ipc.WINAC.Support1InB'  
in_2_1 = 'hvl-ipc.WINAC.Support2InA'  
in_2_2 = 'hvl-ipc.WINAC.Support2InB'  
in_3_1 = 'hvl-ipc.WINAC.Support3InA'  
in_3_2 = 'hvl-ipc.WINAC.Support3InB'  
in_4_1 = 'hvl-ipc.WINAC.Support4InA'  
in_4_2 = 'hvl-ipc.WINAC.Support4InB'  
in_5_1 = 'hvl-ipc.WINAC.Support5InA'  
in_5_2 = 'hvl-ipc.WINAC.Support5InB'  
in_6_1 = 'hvl-ipc.WINAC.Support6InA'  
in_6_2 = 'hvl-ipc.WINAC.Support6InB'  
classmethod input(port, contact)
```

Get the NodeID string for a support input.

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string

```
out_1_1 = 'hvl-ipc.WINAC.Support1OutA'  
out_1_2 = 'hvl-ipc.WINAC.Support1OutB'  
out_2_1 = 'hvl-ipc.WINAC.Support2OutA'  
out_2_2 = 'hvl-ipc.WINAC.Support2OutB'  
out_3_1 = 'hvl-ipc.WINAC.Support3OutA'  
out_3_2 = 'hvl-ipc.WINAC.Support3OutB'  
out_4_1 = 'hvl-ipc.WINAC.Support4OutA'  
out_4_2 = 'hvl-ipc.WINAC.Support4OutB'  
out_5_1 = 'hvl-ipc.WINAC.Support5OutA'  
out_5_2 = 'hvl-ipc.WINAC.Support5OutB'  
out_6_1 = 'hvl-ipc.WINAC.Support6OutA'  
out_6_2 = 'hvl-ipc.WINAC.Support6OutB'
```

```
classmethod output(port, contact)
```

Get the NodeID string for a support output.

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string

```
class MeasurementsDividerRatio(value=<no_arg>, names=None, module=None, type=None, start=1,
                                 boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

classmethod get(channel: int)

Get the attribute for an input number.

Parameters `channel` – the channel number (1..4)

Returns the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.DivididerRatio'
```

```
class MeasurementsScaledInput(value=<no_arg>, names=None, module=None, type=None, start=1,
                                 boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in `MeasurementsDividerRatio`.

classmethod get(channel: int)

Get the attribute for an input number.

Parameters `channel` – the channel number (1..4)

Returns the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI1Volt'
```

```
input_2 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI2Volt'
```

```
input_3 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI3Volt'
```

```
input_4 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI4Volt'
```

```
class Power(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

Variable NodeID strings concerning power data.

```
current_primary = 'hvl-ipc.WINAC.SYSTEM_INTERN.FUCurrentprim'
```

Primary current in ampere, measured by the frequency converter. (read-only)

```
frequency = 'hvl-ipc.WINAC.FU.Frequency'
```

Frequency converter output frequency. (read-only)

```
setup = 'hvl-ipc.WINAC.FU.TrafoSetup'
```

Power setup that is configured using the Supercube HMI. The value corresponds to the ones in `PowerSetup`. (read-only)

```
voltage_max = 'hvl-ipc.WINAC.FU.maxVoltagekV'
```

Maximum voltage allowed by the current experimental setup. (read-only)

```
voltage_primary = 'hvl-ipc.WINAC.SYSTEM_INTERN.FUVoltageprim'
```

Primary voltage in volts, measured by the frequency converter at its output. (read-only)

```
voltage_slope = 'hvl-ipc.WINAC.FU.dUdt_-1'
```

Voltage slope in V/s.

```
voltage_target = 'hvl-ipc.WINAC.FU.SOLL'
```

Target voltage setpoint in V.

```
class PowerSetup(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `aenum.IntEnum`

Possible power setups corresponding to the value of variable `Power.setup`.

AC_DoubleStage_150kV = 3

AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum voltage of 150kV.

AC_DoubleStage_200kV = 4

AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

AC_SingleStage_100kV = 2

AC voltage with MWB transformer set to 100kV maximum voltage.

AC_SingleStage_50kV = 1

AC voltage with MWB transformer set to 50kV maximum voltage.

DC_DoubleStage_280kV = 7

DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single stage transformer with Greinacher voltage doubling rectifier)

DC_SingleStage_140kV = 6

DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

External = 0

External power supply fed through blue CEE32 input using isolation transformer and safety switches of the Supercube, or using an external safety switch attached to the Supercube Type B.

Internal = 5

Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

```
class Safety(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

horn = 'hvl-ipc.WINAC.SYSTEM_INTERN.hornen'

Writeable boolean to manually turn on or off the horn

status_error = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'

status_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN'

status_ready_for_red = 'hvl-ipc.WINAC.SYSTEMSTATE.ReadyForRed'

Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in `SafetyStatus`.

status_red = 'hvl-ipc.WINAC.SYSTEMSTATE.RED'

switchto_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN_REQUEST'

switchto_operate = 'hvl-ipc.WINAC.SYSTEMSTATE.switchon'

Writable boolean for switching to Red Operate (locket, HV on) state.

switchto_ready = 'hvl-ipc.WINAC.SYSTEMSTATE.RED_REQUEST'

Writable boolean for switching to Red Ready (locked, HV off) state.

```
class SafetyStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `aenum.IntEnum`

Safety status values that are possible states returned from `hvl_ccb.dev.supercube.base.Supercube.get_status()`. These values correspond to the states of the Supercube’s safety circuit statemachine.

Error = 6

System is in error mode.

GreenNotReady = 1

System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

GreenReady = 2

System is safe and all safety elements are in place to be able to switch to *ready*.

Initializing = 0

System is initializing or booting.

QuickStop = 5

Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

RedOperate = 4

System is locked in red state and in *operate* mode, i.e. high voltage on.

RedReady = 3

System is locked in red state and *ready* to go to *operate* mode.

```
class SupercubeOpcEndpoint(value=<no_arg>, names=None, module=None, type=None, start=1,  
                             boundary=None)
```

Bases: *hvl_ccb.utils.enum.ValueEnum*

OPC Server Endpoint strings for the supercube variants.


A = 'OPC.SimaticNET.S7'

B = 'OPC.SimaticNET.S7'

```
T13_SOCKET_PORTS = (1, 2, 3)
```

Port numbers of SEV T13 power socket

`hvl_ccb.dev.supercube2015.typ_a`

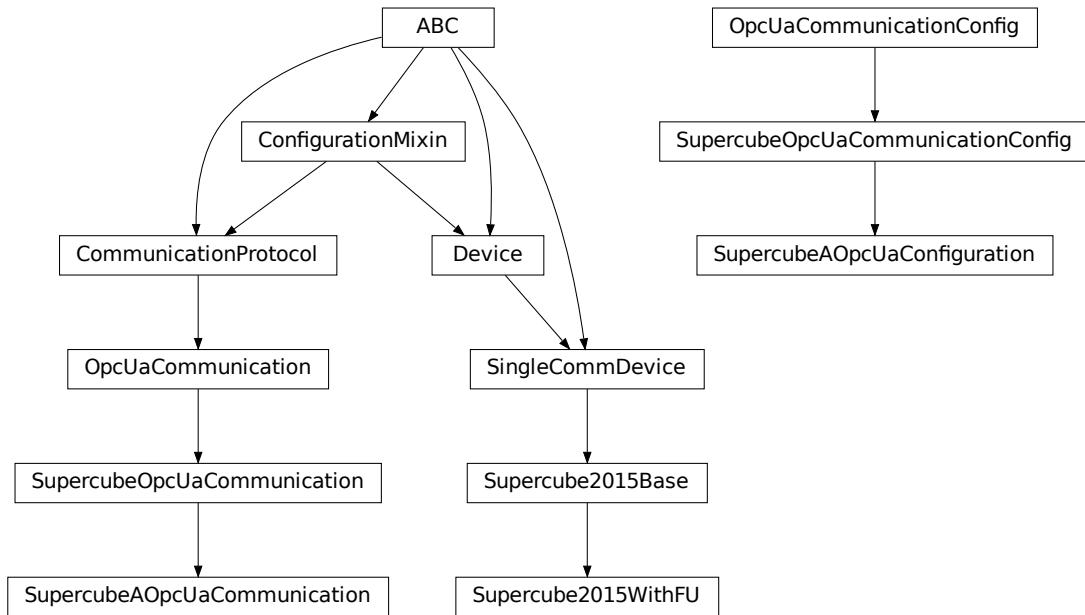

Supercube Typ A module.

class Supercube2015WithFU(com, dev_config=None)
Bases: `hvl_ccb.dev.supercube2015.base.Supercube2015Base`

Variant A of the Supercube with frequency converter.

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

fso_reset() → None

Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.

get_frequency() → float

Read the electrical frequency of the current Supercube setup.

Returns the frequency in Hz

get_fso_active() → bool

Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.

Returns state of the FSO functionality

get_max_voltage() → float

Reads the maximum voltage of the setup and returns in V.

Returns the maximum voltage of the setup in V.

get_power_setup() → `hvl_ccb.dev.supercube2015.constants.PowerSetup`

Return the power setup selected in the Supercube's settings.

Returns the power setup

get_primary_current() → float
Read the current primary current at the output of the frequency converter (before transformer).

Returns primary current in A

get_primary_voltage() → float
Read the current primary voltage at the output of the frequency converter (before transformer).

Returns primary voltage in V

get_target_voltage() → float
Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to `set_target_voltage()`.

Returns the setpoint voltage in V.

set_slope(*slope: float*) → None
Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.

Parameters **slope** – voltage slope in V/s (0..15'000)

set_target_voltage(*volt_v: float*) → None
Set the output voltage to a defined value in V.

Parameters **volt_v** – the desired voltage in V

class SupercubeA0pcUaCommunication(*config*)
Bases: `hvl_ccb.dev.supercube2015.base.SupercubeOpcUaCommunication`

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

class SupercubeA0pcUaConfiguration(*host: str, endpoint_name: str = 'OPC.SimaticNET.S7', port: int = 4845, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object at 0x7fed1b06a4d0>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5*)
Bases: `hvl_ccb.dev.supercube2015.base.SupercubeOpcUaCommunicationConfig`

endpoint_name: str = 'OPC.SimaticNET.S7'
Endpoint of the OPC server, this is a path like ‘OPCUA/SimulationServer’

force_value(*fieldname, value*)
Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.


Module contents

Supercube package with implementation for the old system version from 2015 based on Siemens WinAC soft-PLC on an industrial 32bit Windows computer.

[hvl_ccb.dev.tiepie](#)

Submodules

[hvl_ccb.dev.tiepie.base](#)

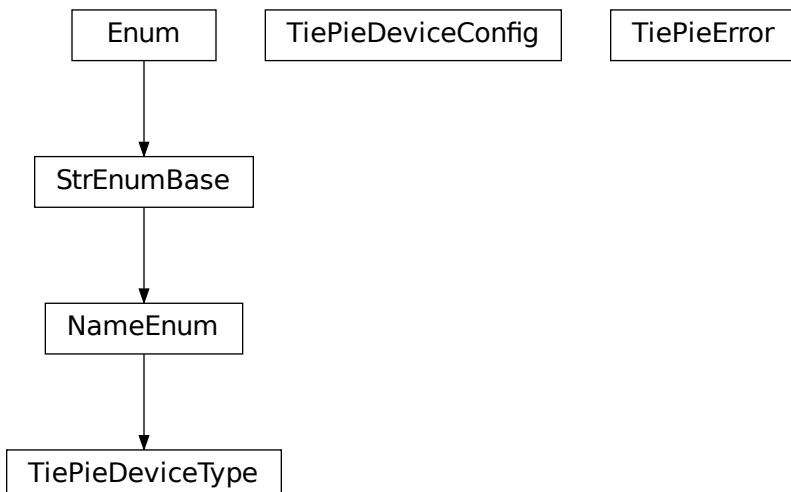

class TiePieDeviceConfig(*serial_number: int, require_block_measurement_support: bool = True, n_max_try_get_device: int = 10, wait_sec_retry_get_device: Union[int, float] = 1.0, is_data_ready_polling_interval_sec: Union[int, float] = 0.01*)

Bases: `object`

Configuration dataclass for TiePie

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**is_data_ready_polling_interval_sec: Union[int, float] = 0.01****classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

n_max_try_get_device: int = 10**classmethod optional_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

require_block_measurement_support: bool = True**classmethod required_keys()** → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

serial_number: int**wait_sec_retry_get_device: Union[int, float] = 1.0****class TiePieDeviceType**(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: *hvl_ccb.utils.enum.NameEnum*

TiePie device type.

GENERATOR = 2**I2C = 4****OSCILLOSCOPE = 1****exception TiePieError**

Bases: *Exception*

Error of the class TiePie

get_device_by_serial_number(*serial_number: int*, *device_type: Union[str, Tuple[int, hvl_ccb.dev.tiepie.base._LtpDeviceReturnType]]*, *n_max_try_get_device: int = 10*, *wait_sec_retry_get_device: float = 1.0*) → *hvl_ccb.dev.tiepie.base._LtpDeviceReturnType*

Open and return handle of TiePie device with a given serial number

Parameters

- **serial_number** – int serial number of the device

- **device_type** – a *TiePieDeviceType* instance containing device identifier (int number) and its corresponding class, both from *libtiepie*, or a string name of such instance
- **n_max_try_get_device** – maximal number of device list updates (int number)
- **wait_sec_retry_get_device** – waiting time in seconds between retries (int number)

Returns Instance of a *libtiepie* device class according to the specified *device_type*

Raises

- **TiePieError** – when there is no device with given serial number
- **ValueError** – when *device_type* is not an instance of *TiePieDeviceType*

wrap_libtiepie_exception(func: Callable) → Callable


Decorator wrapper for *libtiepie* methods that use *libtiepie.library.check_last_status_raise_on_error()* calls.

Parameters **func** – Function or method to be wrapped

Raises **TiePieError** – instead of *LibTiePieException* or one of its subtypes.

Returns whatever *func* returns

hvl_ccb.dev.tiepie.channel

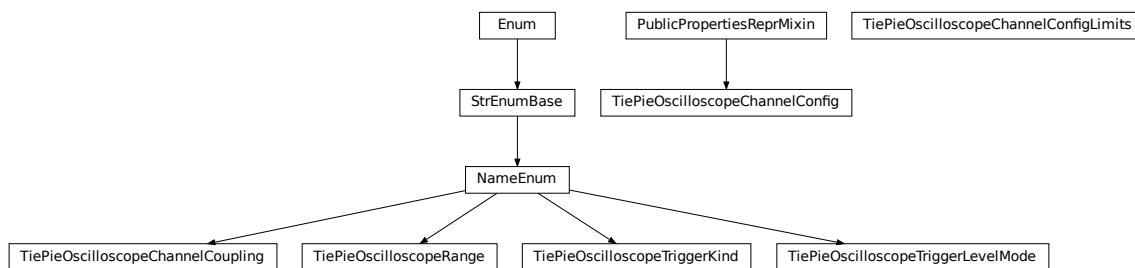

```

class TiePieOscilloscopeChannelConfig(ch_number: int, channel:
libtiepie.oscilloscopechannel.OscilloscopeChannel)
Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin

Oscilloscope's channel configuration, with cleaning of values in properties setters as well as setting and reading them on and from the device's channel.

static clean_coupling(coupling: Union[str,
hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling]) →
hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling

static clean_enabled(enabled: bool) → bool
clean_input_range(input_range: Union[float, hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeRange]) →
hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeRange

clean_probe_offset(probe_offset: float) → float
static clean_safe_ground_enabled(safe_ground_enabled: bool) → bool
static clean_trigger_enabled(trigger_enabled)
  
```

```

clean_trigger_hysteresis(trigger_hysteresis: float) → float
static clean_trigger_kind(trigger_kind: Union[str,
                                              hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind]) →
                                              hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind

clean_trigger_level(trigger_level: Union[int, float]) → float
static clean_trigger_level_mode(level_mode: Union[str,
                                                    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode]) →
                                                    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode

property coupling: hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling
property enabled: bool
property has_safe_ground: bool
    Check whether bound oscilloscope device has “safe ground” option
        Returns bool: 1=safe ground available
property input_range: hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeRange
property probe_offset: float
property safe_ground_enabled: Optional[bool]
property trigger_enabled: bool
property trigger_hysteresis: float
property trigger_kind: hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind
property trigger_level: float
property trigger_level_mode:
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode

class TiePieOscilloscopeChannelConfigLimits(osc_channel:
                                                libtiepie.oscilloscopechannel.OscilloscopeChannel)
Bases: object
Default limits for oscilloscope channel parameters.

class TiePieOscilloscopeChannelCoupling(value=<no_arg>, names=None, module=None, type=None,
                                            start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
An enumeration.

ACA = 8
ACV = 2
DCA = 4
DCV = 1

class TiePieOscilloscopeRange(value=<no_arg>, names=None, module=None, type=None, start=1,
                                 boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
An enumeration.

EIGHTY_VOLT = 80
EIGHT_HUNDRED_MILLI_VOLT = 0.8

```


```
EIGHT_VOLT = 8
FORTY_VOLT = 40
FOUR_HUNDRED_MILLI_VOLT = 0.4
FOUR_VOLT = 4
TWENTY_VOLT = 20
TWO_HUNDRED_MILLI_VOLT = 0.2
TWO_VOLT = 2
static suitable_range(value)

class TiePie0scopeTriggerKind(value=<no_arg>, names=None, module=None, type=None, start=1,
                               boundary=None)
Bases: hvl\_ccb.utils.enum.NameEnum
An enumeration.

ANY = 16
FALLING = 2
RISING = 1
RISING_OR_FALLING = 16

class TiePie0scopeTriggerLevelMode(value=<no_arg>, names=None, module=None, type=None,
                                    start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.NameEnum
An enumeration.

ABSOLUTE = 2
RELATIVE = 1
UNKNOWN = 0
```
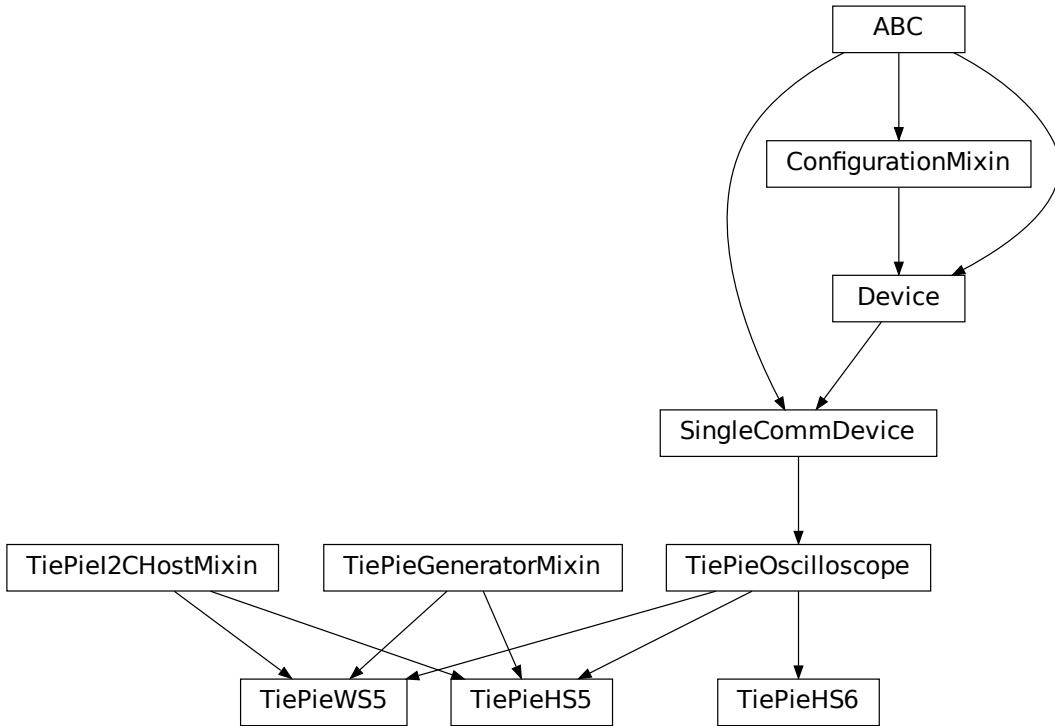
`hvl_ccb.dev.tiepie.device`

TiePie devices.

```
class TiePieHS5(com, dev_config)
Bases:      hvl_ccb.dev.tiepie.i2c.TiePieI2CHostMixin,  hvl_ccb.dev.tiepie.generator.
           TiePieGeneratorMixin, hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope
```

TiePie HS5 device.


```
class TiePieHS6(com, dev_config)
Bases: hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope
```

TiePie HS6 DIFF device.

```
class TiePieWS5(com, dev_config)
Bases:      hvl_ccb.dev.tiepie.i2c.TiePieI2CHostMixin,  hvl_ccb.dev.tiepie.generator.
           TiePieGeneratorMixin, hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope
```

TiePie WS5 device.

`hvl_ccb.dev.tiepie.generator`

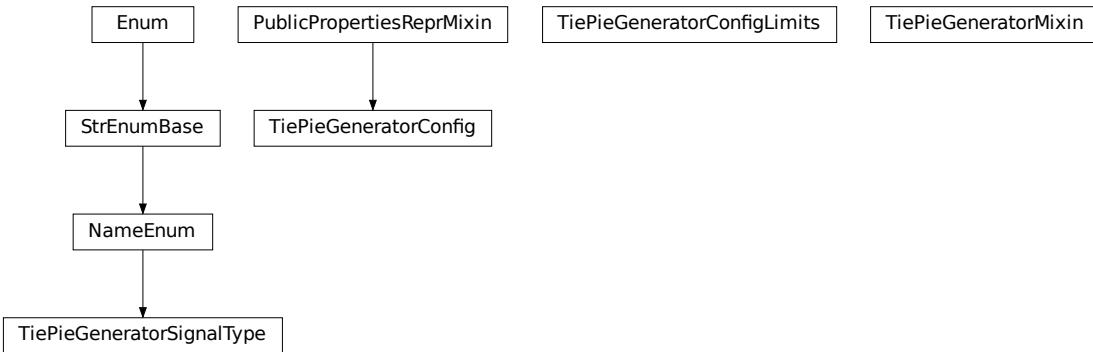

```

class TiePieGeneratorConfig(dev_gen: libtiepie.generator.Generator)
  Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin
  Generator's configuration with cleaning of values in properties setters.

  property amplitude: float
  clean_amplitude(amplitude: float) → float
  static clean_enabled(enabled: bool) → bool
  clean_frequency(frequency: float) → float
  clean_offset(offset: float) → float
  static clean_signal_type(signal_type: Union[int,
                                              hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType]) →
                                              hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType

  property enabled: bool
  property frequency: float
  property offset: float
  property signal_type: hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType

class TiePieGeneratorConfigLimits(dev_gen: libtiepie.generator.Generator)
  Bases: object
  Default limits for generator parameters.

class TiePieGeneratorMixin(com, dev_config)
  Bases: object
  TiePie Generator sub-device.

  A wrapper for the libtiepie.generator.Generator class. To be mixed in with TiePieOscilloscope base class.

  config_gen: Optional[hvl_ccb.dev.tiepie.generator.TiePieGeneratorConfig]
  Generator's dynamical configuration.
  
```

```

generator_start()
    Start signal generation.

generator_stop()
    Stop signal generation.

start() → None
    Start the Generator.

stop() → None
    Stop the generator.

class TiePieGeneratorSignalType(value=<no_arg>, names=None, module=None, type=None, start=1,  


                                    boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ARBITRARY = 32
DC = 8
NOISE = 16
PULSE = 64
SINE = 1
SQUARE = 4
TRIANGLE = 2
UNKNOWN = 0

```

hvl_ccb.dev.tiepie.i2c

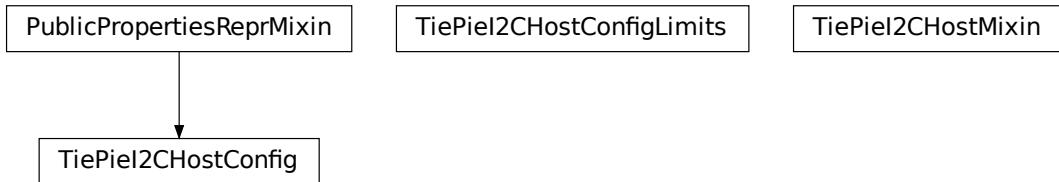

```

class TiePieI2CHostConfig(dev_i2c: libtiepie.i2chost.I2CHost)
Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin

I2C Host's configuration with cleaning of values in properties setters.

class TiePieI2CHostConfigLimits(dev_i2c: libtiepie.i2chost.I2CHost)
Bases: object

Default limits for I2C host parameters.

```

```
class TiePieI2CHostMixin(com, dev_config)
```

Bases: object

TiePie I2CHost sub-device.

A wrapper for the *libtiepie.i2chost.I2CHost* class. To be mixed in with *TiePieOscilloscope* base class.

```
config_i2c: Optional[hvl_ccb.dev.tiepie.i2c.TiePieI2CHostConfig]
```

I2C host's dynamical configuration.


```
start() → None
```

Start the I2C Host.

```
stop() → None
```

Stop the I2C host.

hvl_ccb.dev.tiepie.oscilloscope

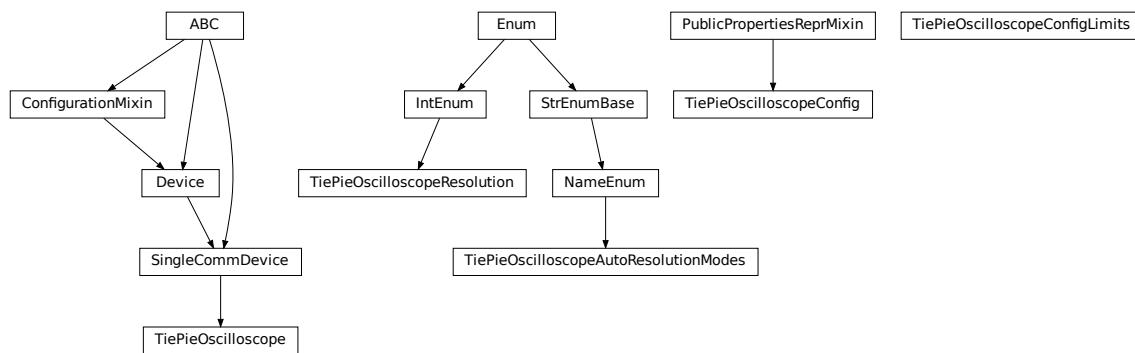

```
class TiePieOscilloscope(com, dev_config)
```

Bases: [hvl_ccb.dev.base.SingleCommDevice](#)

TiePie oscilloscope.

A wrapper for TiePie oscilloscopes, based on the class *libtiepie.oscilloscope.Oscilloscope* with simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and its channels, including extra validation and typing hints support for configurations.

Note that, in contrast to *libtiepie* library, since all physical TiePie devices include an oscilloscope, this is the base class for all physical TiePie devices. The additional TiePie sub-devices: “Generator” and “I2CHost”, are mixed-in to this base class in subclasses.

The channels use *1..N* numbering (not *0..N-1*), as in, e.g., the Multi Channel software.

```
property channels_enabled: Generator[int, None, None]
```

Yield numbers of enabled channels.

Returns Numbers of enabled channels

```
collect_measurement_data(timeout: Optional[Union[int, float]] = 0) → Optional[numpy.ndarray]
```

Try to collect the data from TiePie; return *None* if data is not ready.

Parameters `timeout` – The timeout to wait until data is available. This option makes this function blocking the code. `timeout = None` blocks the code infinitely till data will be available. Per default, the `timeout` is set to `0`: The function will not block.

Returns Measurement data of only enabled channels and time vector in a 2D-`numpy.ndarray` with float sample data; or `None` if there is no data available.

static config_cls() → Type[[hvl_ccb.dev.tiepie.base.TiePieDeviceConfig](#)]
Return the default configdataclass class.

Returns a reference to the default configdataclass class

config_osc: `Optional[hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeConfig]`
Oscilloscope's dynamical configuration.

config_osc_channel_dict: `Dict[int, hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelConfig]`
Channel configuration. A `dict` mapping actual channel number, numbered `1..N`, to channel configuration. The channel info is dynamically read from the device only on the first `start()`; beforehand the `dict` is empty.

static default_com_cls() → Type[[hvl_ccb.comm.base.NullCommunicationProtocol](#)]
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

force_trigger() → `None`
Forces the TiePie to trigger with a software sided trigger event.

Return None

Raises `TiePieError` – when device is not started or status of underlying device gives an error

is_measurement_data_ready() → `bool`
Reports if TiePie has data which is ready to collect

Returns if the data is ready to collect.

Raises `TiePieError` – when device is not started or status of underlying device gives an error

is_triggered() → `bool`
Reports if TiePie has triggered. Maybe data is not yet available. One can check with the function `is_measurement_data_ready()`.

Returns if a trigger event occurred

static list_devices() → libtiepie.devicelist.DeviceList
List available TiePie devices.

Returns libtiepie up to date list of devices

property n_channels
Number of channels in the oscilloscope.

Returns Number of channels.

start() → `None`
Start the oscilloscope.

start_measurement() → `None`
Start a measurement using set configuration.

Raises `TiePieError` – when device is not started or status of underlying device gives an error

stop() → `None`
Stop the oscilloscope.

```
class TiePieOscilloscopeAutoResolutionModes(value=<no_arg>, names=None, module=None,
                                             type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ALL = 4
DISABLED = 1
NATIVEONLY = 2
UNKNOWN = 0

class TiePieOscilloscopeConfig(dev_osc: libtiepie.oscilloscope.Oscilloscope)
Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin

Oscilloscope's configuration with cleaning of values in properties setters.

property auto_resolution_mode:
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes

static clean_auto_resolution_mode(auto_resolution_mode: Union[int,
                                                               hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes])
    →
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes

clean_pre_sample_ratio(pre_sample_ratio: float) → float
clean_record_length(record_length: Union[int, float]) → int
static clean_resolution(resolution: Union[int,
                                            hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution]) →
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution

clean_sample_frequency(sample_frequency: float) → float
clean_trigger_timeout(trigger_timeout: Optional[Union[int, float]]) → float
property pre_sample_ratio: float
property record_length: int
property resolution: hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution
property sample_frequency: float
property trigger_timeout: Optional[float]

class TiePieOscilloscopeConfigLimits(dev_osc: libtiepie.oscilloscope.Oscilloscope)
Bases: object

Default limits for oscilloscope parameters.

class TiePieOscilloscopeResolution(value=<no_arg>, names=None, module=None, type=None, start=1,
                                     boundary=None)
Bases: aenum.IntEnum

An enumeration.

EIGHT_BIT = 8
FOURTEEN_BIT = 14
SIXTEEN_BIT = 16
TWELVE_BIT = 12
```

hvl_ccb.dev.tiepie.utils

PublicPropertiesReprMixin

class PublicPropertiesReprMixin

Bases: object

General purpose utility mixin that overwrites object representation to a one analogous to *dataclass* instances, but using public properties and their values instead of *fields*.

Module contents

This module is a wrapper around LibTiePie SDK devices; see <https://www.tiepie.com/en/libtiepie-sdk> .

The device classes adds simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and oscilloscope's channels. This includes extra validation and typing hints support.

Extra installation

LibTiePie SDK library is available only on Windows and on Linux.

To use this LibTiePie SDK devices wrapper:

1. install the hvl_ccb package with a tiepie extra feature:

```
$ pip install "hvl_ccb[tiepie]"
```

this will install the Python bindings for the library.

2. install the library

- on Linux: follow instructions in <https://www.tiepie.com/en/libtiepie-sdk/linux> ;
- on Windows: the additional DLL is included in Python bindings package.

Troubleshooting

On a Windows system, if you encounter an OSError like this:

```
...
    self._handle = _dlopen(self._name, mode)
OSError: [WinError 126] The specified module could not be found
```


most likely the python-libtiepie package was installed in your site-packages/ directory as a python-libtiepie-*.egg file via python setup.py install or python setup.py develop command. In such case uninstall the library and re-install it using pip:

```
$ pip uninstall python-libtiepie
$ pip install python-libtiepie
```

This should create `libtiepie/` folder. Alternatively, manually move the folder `libtiepie/` from inside of the `.egg` archive file to the containing it `site-packages/` directory (PyCharm's Project tool window supports reading and extracting from `.egg` archives).

Submodules

`hvl_ccb.dev.base`

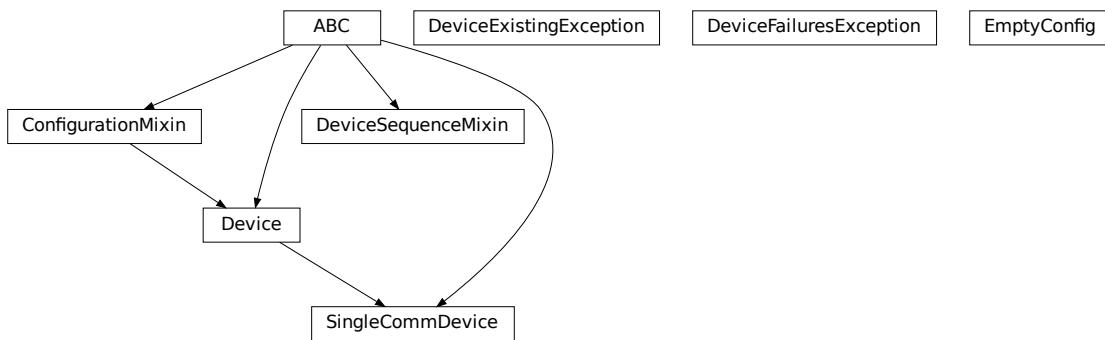

Module with base classes for devices.

class Device(`dev_config=None`)
Bases: `hvl_ccb.configuration.ConfigurationMixin, abc.ABC`

Base class for devices. Implement this class for a concrete device, such as measurement equipment or voltage sources.

Specifies the methods to implement for a device.

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

abstract start() → None
Start or restart this Device. To be implemented in the subclass.

abstract stop() → None
Stop this Device. To be implemented in the subclass.

exception DeviceExistingException
Bases: `Exception`

Exception to indicate that a device with that name already exists.

exception DeviceFailuresException(`failures: Dict[str, Exception], *args`)
Bases: `Exception`

Exception to indicate that one or several devices failed.

failures: Dict[str, Exception]

A dictionary of named devices failures (exceptions).

class DeviceSequenceMixin(devices: Dict[str, hvl_ccb.dev.base.Device])

Bases: abc.ABC

Mixin that can be used on a device or other classes to provide facilities for handling multiple devices in a sequence.

add_device(name: str, device: hvl_ccb.dev.base.Device) → None

Add a new device to the device sequence.

Parameters

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

Raises DeviceExistingException –

devices_failed_start: Dict[str, hvl_ccb.dev.base.Device]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices_failed_stop*.

devices_failed_stop: Dict[str, hvl_ccb.dev.base.Device]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices_failed_start*.

get_device(name: str) → hvl_ccb.dev.base.Device

Get a device by name.

Parameters **name** – is the name of the device.

Returns the device object from this sequence.

get_devices() → List[Tuple[str, hvl_ccb.dev.base.Device]]

Get list of name, device pairs according to current sequence.

Returns A list of tuples with name and device each.

remove_device(name: str) → hvl_ccb.dev.base.Device

Remove a device from this sequence and return the device object.

Parameters **name** – is the name of the device.

Returns device object or *None* if such device was not in the sequence.

Raises ValueError – when device with given name was not found

start() → None

Start all devices in this sequence in their added order.

Raises DeviceFailuresException – if one or several devices failed to start

stop() → None

Stop all devices in this sequence in their reverse order.

Raises DeviceFailuresException – if one or several devices failed to stop

class EmptyConfig

Bases: object

Empty configuration dataclass that is the default configuration for a Device.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

class SingleCommDevice(com, dev_config=None)

Bases: *hvl_ccb.dev.base.Device*, *abc.ABC*

Base class for devices with a single communication protocol.

property com

Get the communication protocol of this device.

Returns an instance of CommunicationProtocol subtype

abstract static default_com_cls() → Type[hvl_ccb.comm.base.CommunicationProtocol]

Get the class for the default communication protocol used with this device.


Returns the type of the standard communication protocol for this device

start() → None

Open the associated communication protocol.

stop() → None

Close the associated communication protocol.

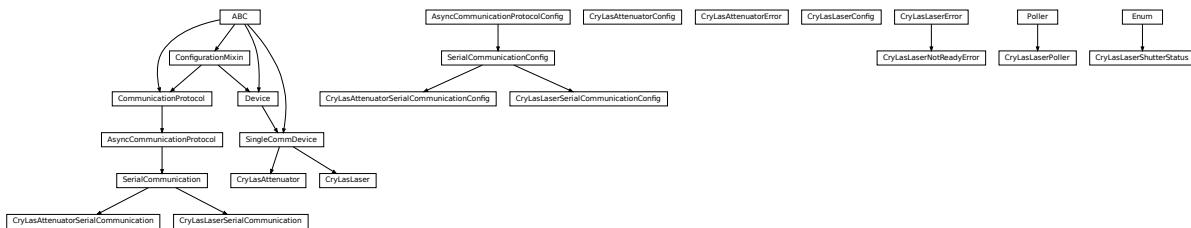hvl_ccb.dev.crylas

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: https://www.crylas.de/products/pulsed_laser.html

class CryLasAttenuator(*com, dev_config=None*)
Bases: *hvl_ccb.dev.base.SingleCommDevice*

Device class for the CryLas laser attenuator.

property attenuation: Union[int, float]

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

set_attenuation(*percent: Union[int, float]*) → None

Set the percentage of attenuated light (inverse of set_transmission). :param percent: percentage of attenuation, number between 0 and 100 :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

set_init_attenuation()

Sets the attenuation to its configured initial/default value

Raises *SerialCommunicationIOError* – when communication port is not opened

set_transmission(*percent: Union[int, float]*) → None

Set the percentage of transmitted light (inverse of set_attenuation). :param percent: percentage of transmitted light :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

start() → None

Open the com, apply the config value ‘init_attenuation’

Raises *SerialCommunicationIOError* – when communication port cannot be opened

property transmission: Union[int, float]

```
class CryLasAttenuatorConfig(init_attenuation: Union[int, float] = 0, response_sleep_time: Union[int, float]
                               = 1)
```

Bases: object

Device configuration dataclass for CryLas attenuator.

clean_values()

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

init_attenuation: Union[int, float] = 0

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

response_sleep_time: Union[int, float] = 1

exception CryLasAttenuatorError

Bases: Exception

General error with the CryLas Attenuator.

class CryLasAttenuatorSerialCommunication(configuration)

Bases: [hvl_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for the CryLas attenuator. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class CryLasAttenuatorSerialCommunicationConfig(terminator: bytes = b'', encoding: str = 'utf-8',
                                                encoding_error_handling: str = 'strict',
                                                wait_sec_read_text_nonempty: Union[int, float] =
                                                0.5, default_n_attempts_read_text_nonempty: int =
                                                10, port: Union[str, NoneType] = None, baudrate: int =
                                                9600, parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for CryLas attenuator is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

CryLas attenuator does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1

CryLas attenuator uses one stop bit

terminator: bytes = b''

No terminator

```
timeout: Union[int, float] = 3
use 3 seconds timeout as default

class CryLasLaser(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice
CryLas laser controller device class.

class AnswersShutter(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: aenum.Enum
Standard answers of the CryLas laser controller to 'Shutter' command passed via com.

CLOSED = 'Shutter inaktiv'
OPENED = 'Shutter aktiv'

class AnswersStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: aenum.Enum
Standard answers of the CryLas laser controller to 'STATUS' command passed via com.

ACTIVE = 'STATUS: Laser active'
HEAD = 'STATUS: Head ok'
INACTIVE = 'STATUS: Laser inactive'
READY = 'STATUS: System ready'
TEC1 = 'STATUS: TEC1 Regulation ok'
TEC2 = 'STATUS: TEC2 Regulation ok'

class LaserStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.Enum
Status of the CryLas laser

READY_ACTIVE = 2
READY_INACTIVE = 1
UNREADY_INACTIVE = 0
property is_inactive
property is_ready

class RepetitionRates(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: aenum.IntEnum
Repetition rates for the internal software trigger in Hz

HARDWARE = 0
SOFTWARE_INTERNAL_SIXTY = 60
SOFTWARE_INTERNAL_TEN = 10
SOFTWARE_INTERNAL_TWENTY = 20

ShutterStatus
alias of hvl_ccb.dev.crylas.CryLasLaserShutterStatus
```

close_shutter() → None

Close the laser shutter.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_pulse_energy_and_rate() → Tuple[int, int]

Use the debug mode, return the measured pulse energy and rate.

Returns (energy in micro joule, rate in Hz)

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not answer the query

laser_off() → None

Turn the laser off.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

laser_on() → None

Turn the laser on.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserNotReadyError* – if the laser is not ready to be turned on
- *CryLasLaserError* – if success is not confirmed by the device

open_shutter() → None

Open the laser shutter.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

set_init_shutter_status() → None

Open or close the shutter, to match the configured shutter_status.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

set_pulse_energy(*energy: int*) → None

Sets the energy of pulses (works only with external hardware trigger). Proceed with small energy steps, or the regulation may fail.

Parameters **energy** – energy in micro joule

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not confirm success

set_repetition_rate(*rate: Union[int, hvl_ccb.dev.crylas.CryLasLaser.RepetitionRates]*) → None

Sets the repetition rate of the internal software trigger.

Parameters **rate** – frequency (Hz) as an integer

Raises

- *ValueError* – if rate is not an accepted value in RepetitionRates Enum
- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

start() → None

Opens the communication protocol and configures the device.

Raises *SerialCommunicationIOError* – when communication port cannot be opened

stop() → None

Stops the device and closes the communication protocol.

Raises

- *SerialCommunicationIOError* – if com port is closed unexpectedly
- *CryLasLaserError* – if laser_off() or close_shutter() fail

property target_pulse_energy

update_laser_status() → None

Update the laser status to *LaserStatus.NOT_READY* or *LaserStatus.INACTIVE* or *LaserStatus.ACTIVE*.

Note: laser never explicitly says that it is not ready (*LaserStatus.NOT_READY*) in response to 'STATUS' command. It only says that it is ready (heated-up and implicitly inactive/off) or active (on). If it's not either of these then the answer is *Answers.HEAD*. Moreover, the only time the laser explicitly says that its status is inactive (*Answers.INACTIVE*) is after issuing a 'LASER OFF' command.

Raises *SerialCommunicationIOError* – when communication port is not opened

update_repetition_rate() → None

Query the laser repetition rate.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

update_shutter_status() → None

Update the shutter status (OPENED or CLOSED)

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

`update_target_pulse_energy()` → None

Query the laser pulse energy.

Raises

- `SerialCommunicationIOError` – when communication port is not opened
- `CryLasLaserError` – if success is not confirmed by the device

`wait_until_ready()` → None

Block execution until the laser is ready

Raises `CryLasLaserError` – if the polling thread stops before the laser is ready

```
class CryLasLaserConfig(calibration_factor: Union[int, float] = 4.35, polling_period: Union[int, float] = 12,
                        polling_timeout: Union[int, float] = 300, auto_laser_on: bool = True,
                        init_shutter_status: Union[int, hvl_ccb.dev.crylas.CryLasLaserShutterStatus] =
                        CryLasLaserShutterStatus.CLOSED)
```

Bases: `object`

Device configuration dataclass for the CryLas laser controller.

ShutterStatus

alias of `hvl_ccb.dev.crylas.CryLasLaserShutterStatus`

`auto_laser_on: bool = True`

`calibration_factor: Union[int, float] = 4.35`

`clean_values()`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- `fieldname` – name of the field
- `value` – value to assign

`init_shutter_status: Union[int, hvl_ccb.dev.crylas.CryLasLaserShutterStatus] = 0`

`is_configdataclass = True`

`classmethod keys()` → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

`classmethod optional_defaults()` → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

`polling_period: Union[int, float] = 12`

`polling_timeout: Union[int, float] = 300`

`classmethod required_keys()` → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception CryLasLaserError

Bases: `Exception`

General error with the CryLas Laser.

exception CryLasLaserNotReadyError

Bases: `hvl_ccb.dev.crylas.CryLasLaserError`

Error when trying to turn on the CryLas Laser before it is ready.

class CryLasLaserPoller(*poll_handler*: Callable, *check_handler*: Callable, *check_laser_status_handler*: Callable, *polling_delay_sec*: Union[int, float] = 0, *polling_interval_sec*: Union[int, float] = 1, *polling_timeout_sec*: Optional[Union[int, float]] = None)

Bases: `hvl_ccb.dev.utils.Poller`

Poller class for polling the laser status until the laser is ready.

Raises

- `CryLasLaserError` – if the timeout is reached before the laser is ready
- `SerialCommunicationIOError` – when communication port is closed.

class CryLasLaserSerialCommunication(*configuration*)

Bases: `hvl_ccb.comm.serial.SerialCommunication`

Specific communication protocol implementation for the CryLas laser controller. Already predefines device-specific protocol parameters in config.

READ_TEXT_SKIP_PREFIXES = ('>', 'MODE:')

Prefixes of lines that are skipped when read from the serial port.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(*cmd*: str, *prefix*: str, *post_cmd*: Optional[str] = None) → str

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns the line in question.

Parameters

- `cmd` – query message to send to the device
- `prefix` – start of the line to look for in the device answer
- `post_cmd` – optional additional command to send after the query

Returns line in question as a string

Raises `SerialCommunicationIOError` – when communication port is not opened

query_all(*cmd*: str, *prefix*: str)

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns a list of successive lines starting with prefix.

Parameters

- `cmd` – query message to send to the device
- `prefix` – start of the line to look for in the device answer

Returns line in question as a string

Raises `SerialCommunicationIOError` – when communication port is not opened

read() → str
 Read first line of text from the serial port that does not start with any of `self.READ_TEXT_SKIP_PREFIXES`.

Returns String read from the serial port; '' if there was nothing to read.

Raises `SerialCommunicationIOError` – when communication port is not opened

```
class CryLasLaserSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Union[str, NoneType] = None, baudrate: int = 19200,
                                             parity: Union[str,
                                             hvl_ccb.comm.serial.SerialCommunicationParity] =
                                             <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                             Union[int,
                                             hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                             <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                             Union[int,
                                             hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                             <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                             Union[int, float] = 10)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 19200

Baudrate for CryLas laser is 19200 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

CryLas laser does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

```

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
    CryLas laser uses one stop bit

terminator: bytes = b'\n'
    The terminator is LF


timeout: Union[int, float] = 10
    use 10 seconds timeout as default (a long timeout is needed!)

class CryLasLaserShutterStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
                                    boundary=None)
Bases: aenum.Enum
Status of the CryLas laser shutter

CLOSED = 0
OPENED = 1

```

hvl_ccb.dev.ea_psi9000

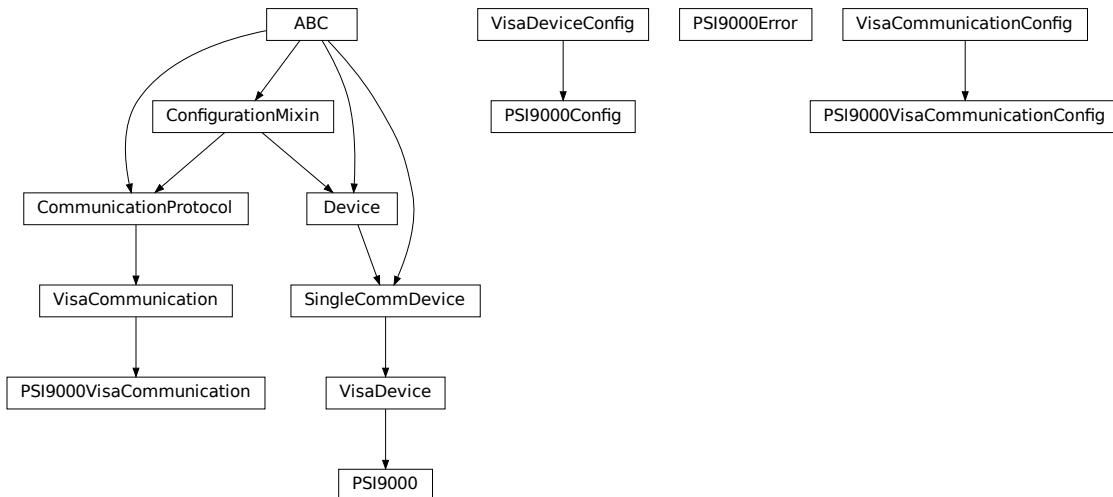

Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to now, all the testing was done with NI-Visa)

```

class PSI9000(com: Union[hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunication,
                    hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunicationConfig, dict], dev_config:
                    Optional[Union[hvl_ccb.dev.ea_psi9000.PSI9000Config, dict]] = None)
Bases: hvl_ccb.dev.visa.VisaDevice
Elektro Automatik PSI 9000 power supply.

MS_NOMINAL_CURRENT = 2040
MS_NOMINAL_VOLTAGE = 80
SHUTDOWN_CURRENT_LIMIT = 0.1

```

SHUTDOWN_VOLTAGE_LIMIT = 0.1

check_master_slave_config() → None
Checks if the master / slave configuration and initializes if successful
Raises *PSI9000Error* – if master-slave configuration failed

static config_cls()
Return the default configdataclass class.
Returns a reference to the default configdataclass class

static default_com_cls()
Return the default communication protocol for this device type, which is VisaCommunication.
Returns the VisaCommunication class

get_output() → bool
Reads the current state of the DC output of the source. Returns True, if it is enabled, false otherwise.
Returns the state of the DC output

get_system_lock() → bool
Get the current lock state of the system. The lock state is true, if the remote control is active and false, if not.
Returns the current lock state of the device

get_ui_lower_limits() → Tuple[float, float]
Get the lower voltage and current limits. A lower power limit does not exist.
Returns Umin in V, Imin in A

get_uip_upper_limits() → Tuple[float, float, float]
Get the upper voltage, current and power limits.
Returns Umax in V, Imax in A, Pmax in W

get_voltage_current_setpoint() → Tuple[float, float]
Get the voltage and current setpoint of the current source.
Returns Uset in V, Iset in A

measure_voltage_current() → Tuple[float, float]
Measure the DC output voltage and current
Returns Umeas in V, Imeas in A

set_lower_limits(voltage_limit: Optional[float] = None, current_limit: Optional[float] = None) → None
Set the lower limits for voltage and current. After writing the values a check is performed if the values are set correctly.

Parameters

- **voltage_limit** – is the lower voltage limit in V
- **current_limit** – is the lower current limit in A

Raises *PSI9000Error* – if the limits are out of range

set_output(target_onstate: bool) → None
Enables / disables the DC output.

Parameters **target_onstate** – enable or disable the output power
Raises *PSI9000Error* – if operation was not successful

set_system_lock(lock: bool) → None

Lock / unlock the device, after locking the control is limited to this class unlocking only possible when voltage and current are below the defined limits

Parameters lock – True: locking, False: unlocking

set_upper_limits(voltage_limit: Optional[float] = None, current_limit: Optional[float] = None, power_limit: Optional[float] = None) → None

Set the upper limits for voltage, current and power. After writing the values a check is performed if the values are set. If a parameter is left blank, the maximum configurable limit is set.

Parameters

- **voltage_limit** – is the voltage limit in V
- **current_limit** – is the current limit in A
- **power_limit** – is the power limit in W

Raises **PSI9000Error** – if limits are out of range

set_voltage_current(volt: float, current: float) → None

Set voltage and current setpoints.

After setting voltage and current, a check is performed if writing was successful.

Parameters

- **volt** – is the setpoint voltage: 0.81.6 V (1.02 * 0-80 V) (absolute max, can be smaller if limits are set)
- **current** – is the setpoint current: 0..2080.8 A (1.02 * 0 - 2040 A) (absolute max, can be smaller if limits are set)

Raises **PSI9000Error** – if the desired setpoint is out of limits

start() → None

Start this device.

stop() → None

Stop this device. Turns off output and lock, if enabled.

class PSI9000Config(spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2, power_limit: Union[int, float] = 43500, voltage_lower_limit: Union[int, float] = 0.0, voltage_upper_limit: Union[int, float] = 10.0, current_lower_limit: Union[int, float] = 0.0, current_upper_limit: Union[int, float] = 2040.0, wait_sec_system_lock: Union[int, float] = 0.5, wait_sec_settings_effect: Union[int, float] = 1, wait_sec_initialisation: Union[int, float] = 2)

Bases: [hvl_ccb.dev.visa.VisaDeviceConfig](#)

Elektro Automatik PSI 9000 power supply device class. The device is communicating over a VISA TCP socket.

Using this power supply, DC voltage and current can be supplied to a load with up to 2040 A and 80 V (using all four available units in parallel). The maximum power is limited by the grid, being at 43.5 kW available through the CEE63 power socket.

clean_values() → None

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

current_lower_limit: Union[int, float] = 0.0

Lower current limit in A, depending on the experimental setup.

current_upper_limit: Union[int, float] = 2040.0

Upper current limit in A, depending on the experimental setup.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

power_limit: Union[int, float] = 43500

Power limit in W depending on the experimental setup. With 3x63A, this is 43.5kW. Do not change this value, if you do not know what you are doing. There is no lower power limit.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

voltage_lower_limit: Union[int, float] = 0.0

Lower voltage limit in V, depending on the experimental setup.

voltage_upper_limit: Union[int, float] = 10.0

Upper voltage limit in V, depending on the experimental setup.

wait_sec_initialisation: Union[int, float] = 2

wait_sec_settings_effect: Union[int, float] = 1

wait_sec_system_lock: Union[int, float] = 0.5

exception PSI9000Error

Bases: Exception

Base error class regarding problems with the PSI 9000 supply.

class PSI9000VisaCommunication(configuration)

Bases: *hvl_ccb.comm.visa.VisaCommunication*

Communication protocol used with the PSI 9000 power supply.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class PSI9000VisaCommunicationConfig(host: str, interface_type: Union[str,
    hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType] =
    InterfaceType.TCPPIP_SOCKET, board: int = 0, port: int = 5025,
    timeout: int = 5000, chunk_size: int = 204800, open_timeout: int =
    1000, write_termination: str = '\n', read_termination: str = '\n',
    visa_backend: str = '')
```

Bases: `hvl_ccb.comm.visa.VisaCommunicationConfig`

Visa communication protocol config dataclass with specification for the PSI 9000 power supply.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

interface_type: `Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]`
= 1

Interface type of the VISA connection, being one of `InterfaceType`.

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.


Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

`hvl_ccb.dev.fug`

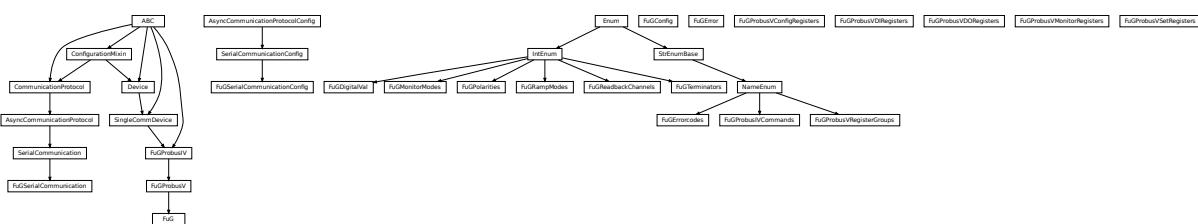

Device classes for “Probus V - ADDAT30” Interfaces which are used to control power supplies from FuG Elektronik GmbH

This interface is used for many FuG power units. Manufacturer homepage: <https://www.fug-elektronik.de>

The Professional Series of Power Supplies from FuG is a series of low, medium and high voltage direct current power supplies as well as capacitor chargers. The class `FuG` is tested with a HCK 800-20 000 in Standard Mode. The

addressable mode is not implemented. Check the code carefully before using it with other devices. Manufacturer homepage: <https://www.fug-elektronik.de/netzgeraete/professional-series/>

The documentation of the interface from the manufacturer can be found here: https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus_V.zip

The provided classes support the basic and some advanced commands. The commands for calibrating the power supplies are not implemented, as they are only for very special purposes and should not be used by “normal” customers.

```
class FuG(com, dev_config=None)
```

Bases: `hvl_ccb.dev.fug.FuGProbusV`

FuG power supply device class.

The power supply is controlled over a FuG ADDA Interface with the PROBUS V protocol

```
property config_status: hvl_ccb.dev.fug.FuGProbusVConfigRegisters
```

Returns the registers for the registers with the configuration and status values

Returns FuGProbusVConfigRegisters

```
property current: hvl_ccb.dev.fug.FuGProbusVSetRegisters
```

Returns the registers for the current output

Returns

```
property current_monitor: hvl_ccb.dev.fug.FuGProbusVMonitorRegisters
```

Returns the registers for the current monitor.

A typical usage will be “self.current_monitor.value” to measure the output current

Returns

```
property di: hvl_ccb.dev.fug.FuGProbusVDIRegisters
```

Returns the registers for the digital inputs

Returns FuGProbusVDIRegisters

```
identify_device() → None
```

Identify the device nominal voltage and current based on its model number.

Raises `SerialCommunicationIOError` – when communication port is not opened

```
property max_current: Union[int, float]
```

Returns the maximal current which could be provided within the test setup

Returns

```
property max_current_hardware: Union[int, float]
```

Returns the maximal current which could be provided with the power supply

Returns

```
property max_voltage: Union[int, float]
```

Returns the maximal voltage which could be provided within the test setup

Returns

```
property max_voltage_hardware: Union[int, float]
```

Returns the maximal voltage which could be provided with the power supply

Returns

```
property on: hvl_ccb.dev.fug.FuGProbusVDORegisters
```

Returns the registers for the output switch to turn the output on or off

Returns FuGProbusVDORegisters

property outX0: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X0

Returns FuGProbusVDORegisters

property outX1: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X1

Returns FuGProbusVDORegisters

property outX2: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X2

Returns FuGProbusVDORegisters

property outXCMD: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital outputX-CMD

Returns FuGProbusVDORegisters

start(*max_voltage=0, max_current=0*) → None

Opens the communication protocol and configures the device.

Parameters

- **max_voltage** – Configure here the maximal permissible voltage which is allowed in the given experimental setup
- **max_current** – Configure here the maximal permissible current which is allowed in the given experimental setup

property voltage: *hvl_ccb.dev.fug.FuGProbusVSetRegisters*

Returns the registers for the voltage output

Returns

property voltage_monitor: *hvl_ccb.dev.fug.FuGProbusVMonitorRegisters*

Returns the registers for the voltage monitor.

A typically usage will be “self.voltage_monitor.value” to measure the output voltage

Returns

class FuGConfig(*wait_sec_stop_commands: Union[int, float] = 0.5*)

Bases: object

Device configuration dataclass for FuG power supplies.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
 Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
 Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

wait_sec_stop_commands: Union[int, float] = 0.5
 Time to wait after subsequent commands during stop (in seconds)

class FuGDigitalVal(value)
 Bases: `enum.IntEnum`

An enumeration.

NO = 0

OFF = 0

ON = 1

YES = 1

exception FuGError(*args, **kwargs)
 Bases: `Exception`

Error with the FuG voltage source.

errorcode: str
 Errorcode from the Probus, see documentation of Probus V chapter 5. Errors with three-digit errorcodes are thrown by this python module.

class FuGErrorcodes(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
 Bases: `hvl_ccb.utils.enum.NameEnum`

The power supply can return an errorcode. These errorcodes are handled by this class. The original errorcodes from the source are with one or two digits, see documentation of Probus V chapter 5. All three-digit errorcodes are from this python module.

E0 = ('no error', 'standard response on each command')

E1 = ('no data available', 'Customer tried to read from GPIB but there were no data prepared. (IBIG50 sent command ~T2 to ADDA)')

E10 = ('unknown SCPI command', 'This SCPI command is not implemented')

E100 = ('Command is not implemented', 'You tried to execute a command, which is not implemented or does not exist')

E106 = ('The rampstate is a read-only register', 'You tried to write data to the register, which can only give you the status of the ramping.')

E11 = ('not allowed Trigger-on-Talk', 'Not allowed attempt to Trigger-on-Talk (~T1) while ADDA was in addressable mode.')

```
E115 = ('The given index to select a digital value is out of range', 'Only integer values between 0 and 1 are allowed.')
E12 = ('invalid argument in ~Tn command', 'Only ~T1 and ~T2 is implemented.')
E125 = ('The given index to select a ramp mode is out of range', 'Only integer values between 0 and 4 are allowed.')
E13 = ('invalid N-value', 'Register > K8 contained an invalid value. Error code is output on an attempt to query data with ? or ~T1')
E135 = ('The given index to select the readback channel is out of range', 'Only integer values between 0 and 6 are allowed.')
E14 = ('register is write only', 'Some registers can only be written to (i.e. > H0)')
E145 = ('The given value for the AD-conversion is unknown', 'Valid values for the ad-conversion are integer values from "0" to "7".')
E15 = ('string too long', 'i.e. serial number string too long during calibration')
E155 = ('The given value to select a polarity is out range.', 'The value should be 0 or 1.')
E16 = ('wrong checksum', 'checksum over command string was not correct, refer also to 4.4 of the Probus V documentation')
E165 = ('The given index to select the terminator string is out of range', '')
E2 = ('unknown register type', "No valid register type after '>'")
E206 = ('This status register is read-only', 'You tried to write data to this register, which can only give you the actual status of the corresponding digital output.')
E306 = ('The monitor register is read-only', 'You tried to write data to a monitor, which can only give you measured data.')
E4 = ('invalid argument', 'The argument of the command was rejected .i.e. malformed number')
E5 = ('argument out of range', 'i.e. setvalue higher than type value')
E504 = ('Empty string as response', 'The connection is broken.')
E505 = ('The returned register is not the requested.', 'Maybe the connection is overburden.')
E6 = ('register is read only', 'Some registers can only be read but not written to. (i.e. monitor registers)')
E666 = ('You cannot overwrite the most recent error in the interface of the power supply. But, well: You created an error anyway...', '')
E7 = ('Receive Overflow', 'Command string was longer than 50 characters.')
E8 = ('EEPROM is write protected', 'Write attempt to calibration data while the write protection switch was set to write protected.')
E9 = ('address error', 'A non addressed command was sent to ADDA while it was in addressable mode (and vice versa).')
raise_()
```

class FuGMonitorModes(*value*)Bases: `enum.IntEnum`

An enumeration.

T1MS = 1

15 bit + sign, 1 ms integration time

T200MS = 6

typ. 19 bit + sign, 200 ms integration time

T20MS = 3

17 bit + sign, 20 ms integration time

T256US = 0

14 bit + sign, 256 us integration time

T40MS = 4

17 bit + sign, 40 ms integration time

T4MS = 2

15 bit + sign, 4 ms integration time

T800MS = 7

typ. 20 bit + sign, 800 ms integration time

T80MS = 5

typ. 18 bit + sign, 80 ms integration time

class FuGPolarities(*value*)Bases: `enum.IntEnum`

An enumeration.

NEGATIVE = 1**POSITIVE = 0****class FuGProbusIV(*com, dev_config=None*)**Bases: `hvl_ccb.dev.base.SingleCommDevice, abc.ABC`**FuG Probus IV device class****Returns** Sends basic SCPI commands and reads the answer. Only the special commands and PROBUS IV instruction set is implemented.**command**(*command: hvl_ccb.dev.fug.FuGProbusIVCommands, value=None*) → str**Parameters**

- **command** – one of the commands given within `FuGProbusIVCommands`
- **value** – an optional value, depending on the command

Returns a String if a query was performed**static config_cls()**

Return the default configdataclass class.

Returns a reference to the default configdataclass class**static default_com_cls()**

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

output_off() → None
Switch DC voltage output off.

reset() → None
Reset of the interface: All setvalues are set to zero

abstract start()
Open the associated communication protocol.

stop() → None
Close the associated communication protocol.

class FuGProbusIVCommands(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)
Bases: *hvl_ccb.utils.enum.NameEnum*

An enumeration.

ADMODE = ('S', (<enum 'FuGMonitorModes'>, <class 'int'>))

CURRENT = ('I', (<class 'int'>, <class 'float'>))

EXECUTE = ('X', None)

EXECUTEONX = ('G', (<enum 'FuGDigitalVal'>, <class 'int'>))
Wait for "X" to execute pending commands

ID = ('*IDN?', None)

OUTPUT = ('F', (<enum 'FuGDigitalVal'>, <class 'int'>))

POLARITY = ('P', (<enum 'FuGPolarities'>, <class 'int'>))

QUERY = ('?', None)

READBACKCHANNEL = ('N', (<enum 'FuGReadbackChannels'>, <class 'int'>))

RESET = ('=', None)

TERMINATOR = ('Y', (<enum 'FuGTerminators'>, <class 'int'>))

VOLTAGE = ('U', (<class 'int'>, <class 'float'>))

XOUTPUTS = ('R', <class 'int'>)
TODO: the possible values are limited to 0..13

class FuGProbusV(*com*, *dev_config*=None)

Bases: *hvl_ccb.dev.fug.FuGProbusIV*

FuG Probus V class which uses register based commands to control the power supplies

get_register(*register*: str) → str
get the value from a register

Parameters **register** – the register from which the value is requested

Returns the value of the register as a String

set_register(*register*: str, *value*: Union[int, float, str]) → None
generic method to set value to register

Parameters

- **register** – the name of the register to set the value
- **value** – which should be written to the register

```
class FuGProbusVConfigRegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Configuration and Status values, acc. 4.2.5

property execute_on_x: hvl_ccb.dev.fug.FuGDigitalVal
    status of Execute-on-X

        Returns FuGDigitalVal of the status

property most_recent_error: hvl_ccb.dev.fug.FuGErrorcodes
    Reads the Error-Code of the most recent command

        Return FuGError

        Raises FuGError – if code is not “E0”

property readback_data: hvl_ccb.dev.fug.FuGReadbackChannels
    Preselection of readout data for Trigger-on-Talk

        Returns index for the readback channel

property srq_mask: int
    SRQ-Mask, Service-Request Enable status bits for SRQ 0: no SRQ Bit 2: SRQ on change of status to CC
    Bit 1: SRQ on change to CV

        Returns representative integer value

property srq_status: str
    SRQ-Statusbyte output as a decimal number: Bit 2: PS is in CC mode Bit 1: PS is in CV mode

        Returns representative string

property status: str
    Statusbyte as a string of 0/1. Combined status (compatibel to Probus IV), MSB first: Bit 7: I-REG Bit 6:
    V-REG Bit 5: ON-Status Bit 4: 3-Reg Bit 3: X-Stat (polarity) Bit 2: Cal-Mode Bit 1: unused Bit 0: SEL-D

        Returns string of 0/1

property terminator: hvl_ccb.dev.fug.FuGTerminators
    Terminator character for answer strings from ADDA

        Returns FuGTerminators

class FuGProbusVDIRegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Digital Inputs acc. 4.2.4

property analog_control: hvl_ccb.dev.fug.FuGDigitalVal

        Returns shows 1 if power supply is controlled by the analog interface

property calibration_mode: hvl_ccb.dev.fug.FuGDigitalVal

        Returns shows 1 if power supply is in calibration mode

property cc_mode: hvl_ccb.dev.fug.FuGDigitalVal

        Returns shows 1 if power supply is in CC mode

property cv_mode: hvl_ccb.dev.fug.FuGDigitalVal

        Returns shows 1 if power supply is in CV mode

property digital_control: hvl_ccb.dev.fug.FuGDigitalVal

        Returns shows 1 if power supply is digitally controlled
```

```
property on: hvl_ccb.dev.fug.FuGDigitalVal
    Returns shows 1 if power supply ON

property reg_3: hvl_ccb.dev.fug.FuGDigitalVal
    For special applications.

    Returns input from bit 3-REG

property x_stat: hvl_ccb.dev.fug.FuGPolarities
    Returns polarity of HVPS with polarity reversal

class FuGProbusVDORegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Digital outputs acc. 4.2.2

property out: Union[int, hvl_ccb.dev.fug.FuGDigitalVal]
    Status of the output according to the last setting. This can differ from the actual state if output should only
    pulse.

    Returns FuGDigitalVal

property status: hvl_ccb.dev.fug.FuGDigitalVal
    Returns the actual value of output. This can differ from the set value if pulse function is used.

    Returns FuGDigitalVal

class FuGProbusVMonitorRegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Analog monitors acc. 4.2.3

property adc_mode: hvl_ccb.dev.fug.FuGMonitorModes
    The programmed resolution and integration time of the AD converter

    Returns FuGMonitorModes

property value: float
    Value from the monitor.

    Returns a float value in V or A

property value_raw: float
    uncalibrated raw value from AD converter

    Returns float value from ADC

class FuGProbusVRegisterGroups(value=<no_arg>, names=None, module=None, type=None, start=1,
                                 boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

CONFIG = 'K'
INPUT = 'D'
MONITOR_I = 'M1'
MONITOR_V = 'M0'
OUTPUTONCMD = 'BON'
OUTPUTX0 = 'B0'
OUTPUTX1 = 'B1'
```

```

OUTPUTX2 = 'B2'
OUTPUTXCMD = 'BX'
SETCURRENT = 'S1'
SETVOLTAGE = 'S0'

class FuGProbusVSetRegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Setvalue control acc. 4.2.1 for the voltage and the current output

property actualsetvalue: float
    The actual valid set value, which depends on the ramp function.

    Returns actual valid set value

property high_resolution: hvl_ccb.dev.fug.FuGDigitalVal
    Status of the high resolution mode of the output.

        Return 0 normal operation
        Return 1 High Res. Mode

property rampmode: hvl_ccb.dev.fug.FuGRampModes
    The set ramp mode to control the setvalue.

    Returns the mode of the ramp as instance of FuGRampModes

property ramprate: float
    The set ramp rate in V/s.

    Returns ramp rate in V/s

property rampstate: hvl_ccb.dev.fug.FuGDigitalVal
    Status of ramp function.

        Return 0 if final setvalue is reached
        Return 1 if still ramping up

property setvalue: float
    For the voltage or current output this setvalue was programmed.

    Returns the programmed setvalue

class FuGRampModes(value)
Bases: enum.IntEnum

An enumeration.

FOLLOWRAMP = 1
    Follow the ramp up- and downwards

IMMEDIATELY = 0
    Standard mode: no ramp

ONLYUPWARDSOFFTOZERO = 4
    Follow the ramp up- and downwards, if output is OFF set value is zero

RAMPUPWARDS = 2
    Follow the ramp only upwards, downwards immediately

SPECIALRAMPUPWARDS = 3
    Follow a special ramp function only upwards

```

```
class FuGReadbackChannels(value)
```

Bases: `enum.IntEnum`

An enumeration.

```
CURRENT = 1
```

```
FIRMWARE = 5
```

```
RATEDCURRENT = 4
```

```
RATEDVOLTAGE = 3
```

```
SN = 6
```

```
STATUSBYTE = 2
```

```
VOLTAGE = 0
```

```
class FuGSerialCommunication(configuration)
```

Bases: `hvl_ccb.comm.serial.SerialCommunication`

Specific communication protocol implementation for FuG power supplies. Already defines device-specific protocol parameters in config.

```
static config_cls()
```

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
query(command: str) → str
```

Send a command to the interface and handle the status message. Eventually raises an exception.

Parameters `command` – Command to send

Raises `FuGError` – if the connection is broken or the error from the power source itself

Returns Answer from the interface or empty string

```
class FuGSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',  
                                    encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:  
                                    Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int =  
                                    10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity:  
                                    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =  
                                    <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,  
                                    hvl_ccb.comm.serial.SerialCommunicationStopbits] =  
                                    <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,  
                                    hvl_ccb.comm.serial.SerialCommunicationBytesize] =  
                                    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int,  
                                    float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

```
baudrate: int = 9600
```

Baudrate for FuG power supplies is 9600 baud

```
bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
```

One byte is eight bits long

```
default_n_attempts_read_text_nonempty: int = 10
```

default number of attempts to read a non-empty text

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

FuG does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1

FuG uses one stop bit

terminator: bytes = b'\n'

The terminator is LF

timeout: Union[int, float] = 3

use 3 seconds timeout as default

wait_sec_read_text_nonempty: Union[int, float] = 0.5

default time to wait between attempts of reading a non-empty text

class FuGTerminators(value)

Bases: enum.IntEnum

An enumeration.


CR = 3

CRLF = 0

LF = 2

LFCR = 1

hvl_ccb.dev.heinzinger

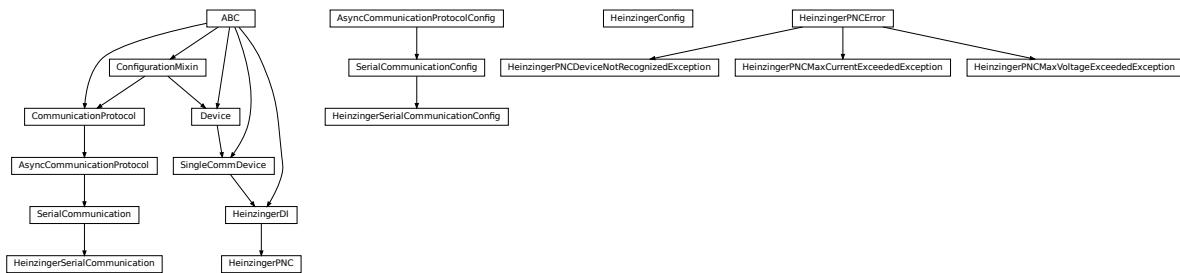

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Manufacturer homepage: <https://www.heinzinger.com/products/accessories-and-more/digital-interfaces/>

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class HeinzingerPNC is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCCap. Manufacturer homepage: <https://www.heinzinger.com/products/high-voltage/universal-high-voltage-power-supplies/>

```

class HeinzingerConfig(default_number_of_recordings: Union[int,
    hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum] = 1,
    number_of_decimals: int = 6, wait_sec_stop_commands: Union[int, float] = 0.5)
  
```

Bases: object

Device configuration dataclass for Heinzinger power supplies.

```

class RecordingsEnum(value)
  
```

Bases: enum. IntEnum

An enumeration.

EIGHT = 8

FOUR = 4

ONE = 1

SIXTEEN = 16

TWO = 2

clean_values()

```

default_number_of_recordings: Union[int,
    hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum] = 1
  
```

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
 Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

number_of_decimals: int = 6

classmethod optional_defaults() → Dict[str, object]
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
 Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

wait_sec_stop_commands: Union[int, float] = 0.5
 Time to wait after subsequent commands during stop (in seconds)

class HeinzingerDI(com, dev_config=None)
 Bases: `hvl_ccb.dev.base.SingleCommDevice`, `abc.ABC`

Heinzinger Digital Interface I/II device class

Sends basic SCPI commands and reads the answer. Only the standard instruction set from the manual is implemented.

class OutputStatus(value)
 Bases: `enum.IntEnum`

Status of the voltage output

OFF = 0

ON = 1

UNKNOWN = -1

static config_cls()
 Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
 Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_current() → float
 Queries the set current of the Heinzinger PNC (not the measured current!).

Raises `SerialCommunicationIOError` – when communication port is not opened

get_interface_version() → str
 Queries the version number of the digital interface.

Raises `SerialCommunicationIOError` – when communication port is not opened

get_number_of_recordings() → int
 Queries the number of recordings the device is using for average value calculation.

Returns int number of recordings

Raises `SerialCommunicationIOError` – when communication port is not opened

`get_serial_number()` → str

Ask the device for its serial number and returns the answer as a string.

Returns string containing the device serial number

Raises `SerialCommunicationIOError` – when communication port is not opened

`get_voltage()` → float

Queries the set voltage of the Heinzinger PNC (not the measured voltage!).

Raises `SerialCommunicationIOError` – when communication port is not opened

`measure_current()` → float

Ask the Device to measure its output current and return the measurement result.

Returns measured current as float

Raises `SerialCommunicationIOError` – when communication port is not opened

`measure_voltage()` → float

Ask the Device to measure its output voltage and return the measurement result.

Returns measured voltage as float

Raises `SerialCommunicationIOError` – when communication port is not opened

`output_off()` → None

Switch DC voltage output off and updates the output status.

Raises `SerialCommunicationIOError` – when communication port is not opened

`output_on()` → None

Switch DC voltage output on and updates the output status.

Raises `SerialCommunicationIOError` – when communication port is not opened

property `output_status: hvl_ccb.dev.heinzinger.HeinzingerDI.OutputStatus`

`reset_interface()` → None

Reset of the digital interface; only Digital Interface I: Power supply is switched to the Local-Mode (Manual operation)

Raises `SerialCommunicationIOError` – when communication port is not opened

`set_current(value: Union[int, float])` → None

Sets the output current of the Heinzinger PNC to the given value.

Parameters `value` – current expressed in `self.unit_current`

Raises `SerialCommunicationIOError` – when communication port is not opened

`set_number_of_recordings(value: Union[int,`

`hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum])` → None

Sets the number of recordings the device is using for average value calculation. The possible values are 1, 2, 4, 8 and 16.

Raises `SerialCommunicationIOError` – when communication port is not opened

`set_voltage(value: Union[int, float])` → None

Sets the output voltage of the Heinzinger PNC to the given value.

Parameters `value` – voltage expressed in `self.unit_voltage`

Raises `SerialCommunicationIOError` – when communication port is not opened

`abstract start()`

Opens the communication protocol.

Raises *SerialCommunicationIOError* – when communication port cannot be opened.

stop() → None

Stop the device. Closes also the communication protocol.

class HeinzingerPNC(*com, dev_config=None*)

Bases: *hvl_ccb.dev.heinzinger.HeinzingerDI*

Heinzinger PNC power supply device class.

The power supply is controlled over a Heinzinger Digital Interface I/II

class UnitCurrent(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl_ccb.utils.enum.AutoNumberNameEnum*

An enumeration.

A = 3

UNKNOWN = 1

mA = 2

class UnitVoltage(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl_ccb.utils.enum.AutoNumberNameEnum*

An enumeration.

UNKNOWN = 1

V = 2

kV = 3

identify_device() → None

Identify the device nominal voltage and current based on its serial number.

Raises *SerialCommunicationIOError* – when communication port is not opened

property max_current: Union[int, float]

property max_current_hardware: Union[int, float]

property max_voltage: Union[int, float]

property max_voltage_hardware: Union[int, float]

set_current(*value: Union[int, float]*) → None

Sets the output current of the Heinzinger PNC to the given value.

Parameters value – current expressed in *self.unit_current*

Raises *SerialCommunicationIOError* – when communication port is not opened

set_voltage(*value: Union[int, float]*) → None

Sets the output voltage of the Heinzinger PNC to the given value.

Parameters value – voltage expressed in *self.unit_voltage*

Raises *SerialCommunicationIOError* – when communication port is not opened

start() → None

Opens the communication protocol and configures the device.

property unit_current: hvl_ccb.dev.heinzinger.HeinzingerPNC.UnitCurrent

property unit_voltage: hvl_ccb.dev.heinzinger.HeinzingerPNC.UnitVoltage

```
exception HeinzingerPNCDeviceNotRecognizedException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError
```

Error indicating that the serial number of the device is not recognized.

```
exception HeinzingerPNCError
```

Bases: Exception

General error with the Heinzinger PNC voltage source.

```
exception HeinzingerPNCMaxCurrentExceededException
```

Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

Error indicating that program attempted to set the current to a value exceeding ‘max_current’.

```
exception HeinzingerPNCMaxVoltageExceededException
```

Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

Error indicating that program attempted to set the voltage to a value exceeding ‘max_voltage’.

```
class HeinzingerSerialCommunication(configuration)
```

Bases: hvl_ccb.comm.serial.SerialCommunication

Specific communication protocol implementation for Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

```
static config_cls()
```

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class HeinzingerSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                            encoding_error_handling: str = 'strict',
                                            wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                            default_n_attempts_read_text_nonempty: int = 40, port:
                                            Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                            Union[str, hvl_ccb.comm.serial.SerialCommunicationParity]
                                            = <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                            Union[int,
                                            hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                            <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                            Union[int,
                                            hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                            <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                            Union[int, float] = 3)
```

Bases: hvl_ccb.comm.serial.SerialCommunicationConfig

```
baudrate: int = 9600
```

Baudrate for Heinzinger power supplies is 9600 baud

```
bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
```

One byte is eight bits long

```
default_n_attempts_read_text_nonempty: int = 40
```

increased to 40 default number of attempts to read a non-empty text

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, *hvl_ccb.comm.serial.SerialCommunicationParity*] = 'N'

Heinzinger does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, *hvl_ccb.comm.serial.SerialCommunicationStopbits*] = 1

Heinzinger uses one stop bit

terminator: bytes = b'\n'

The terminator is LF


timeout: Union[int, float] = 3

use 3 seconds timeout as default

wait_sec_read_text_nonempty: Union[int, float] = 0.5

default time to wait between attempts of reading a non-empty text

hvl_ccb.dev.labjack

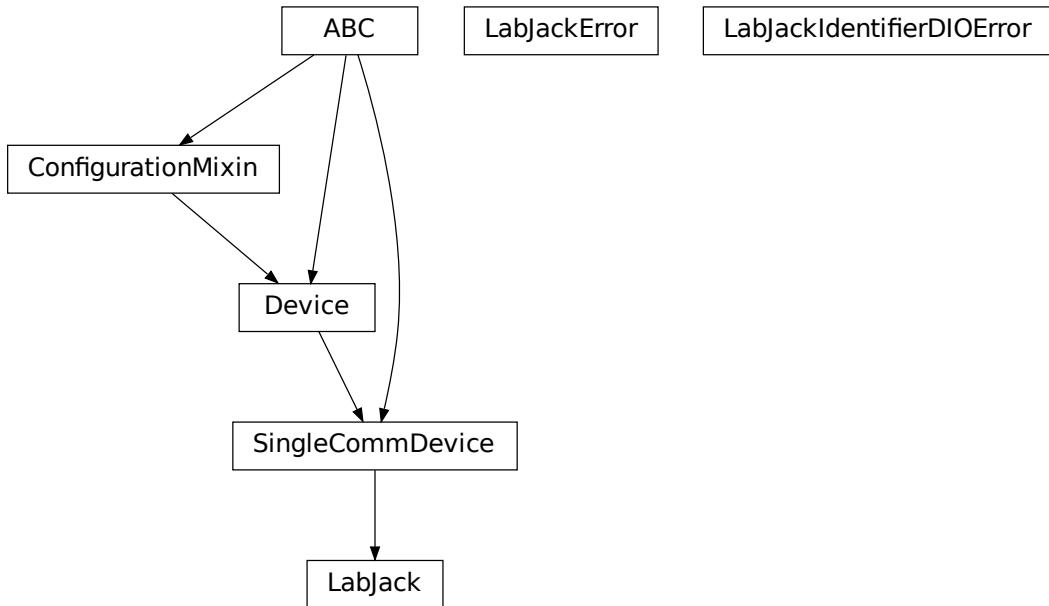

A LabJack T-series devices wrapper around the LabJack's LJM Library; see <https://labjack.com/ljm> . The wrapper was originally developed and tested for a LabJack T7-PRO device.

Extra installation

To use this LabJack T-series devices wrapper:

1. install the hvl_ccb package with a labjack extra feature:

```
$ pip install "hvl_ccb[labjack]"
```

this will install the Python bindings for the library.

2. install the library - follow instruction in <https://labjack.com/support/software/installers/ljm> .

class LabJack(com, dev_config=None)
Bases: *hvl_ccb.dev.base.SingleCommDevice*

LabJack Device.

This class is tested with a LabJack T7-Pro and should also work with T4 and T7 devices communicating through the LJM Library. Other or older hardware versions and variants of LabJack devices are not supported.

class AInRange(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: *hvl_ccb.utils.enum.StrEnumBase*

An enumeration.

ONE = 1.0

```

ONE_HUNDREDTH = 0.01
ONE_TENTH = 0.1
TEN = 10.0
property value: float

class BitLimit(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: aenum.IntEnum
    Maximum integer values for clock settings
    THIRTY_TWO_BIT = 4294967295

class CalMicroAmpere(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
    Bases: aenum.Enum
    Pre-defined microampere (uA) values for calibration current source query.
    TEN = '10uA'
    TWO_HUNDRED = '200uA'

class CjcType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl\_ccb.utils.enum.NameEnum
    CJC slope and offset
    internal = (1, 0)
    lm34 = (55.56, 255.37)

class ClockFrequency(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
    Bases: aenum.IntEnum
    Available clock frequencies, in Hz
    FIVE_MHZ = 5000000
    FORTY_MHZ = 40000000
    MAXIMUM = 80000000
    MINIMUM = 312500
    TEN_MHZ = 10000000
    TWELVE_HUNDRED_FIFTY_KHZ = 1250000
    TWENTY_FIVE_HUNDRED_KHZ = 2500000
    TWENTY_MHZ = 20000000

DIOChannel
    alias of hvl\_ccb.\_dev.labjack.TSeriesDIOChannel

class DIOStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: aenum.IntEnum
    State of a digital I/O channel.
    HIGH = 1
    LOW = 0

```

```
class DeviceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.AutoNumberNameEnum

LabJack device types.

Can be also looked up by ambiguous Product ID (p_id) or by instance name: `python
LabJackDeviceType(4) is LabJackDeviceType('T4')`  

ANY = 1
T4 = 2
T7 = 3
T7_PRO = 4

@classmethod get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
List[hvl_ccb._dev.labjack.DeviceType]]
Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.
Parameters p_id – Product ID of a LabJack device
Returns Instance or list of instances of LabJackDeviceType
Raises ValueError – when Product ID is unknown

class TemperatureUnit(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

Temperature unit (to be returned)

C = 1
F = 2
K = 0

class ThermocoupleType(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

Thermocouple type; NONE means disable thermocouple mode.

C = 30
E = 20
J = 21
K = 22
NONE = 0
PT100 = 40
PT1000 = 42
PT500 = 41
R = 23
S = 25
T = 24
```

config_high_pulse(*address*: Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel], *t_start*: Union[int, float], *t_width*: Union[int, float], *n_pulses*: int = 1) → None

Configures one FIO channel to send a timed HIGH pulse. Configure multiple channels to send pulses with relative timing accuracy. Times have a maximum resolution of 1e-7 seconds @ 10 MHz. :param address: FIO channel: [T7] FIO0;2;3;4;5. [T4] FIO6;7. :raises LabJackError if address is not supported. :param t_start: pulse start time in seconds. :raises ValueError: if t_start is negative or would exceed the clock period. :param t_width: duration of high pulse, in seconds. :raises ValueError: if t_width is negative or would exceed the clock period. :param n_pulses: number of pulses to be sent; single pulse default. :raises TypeError if n_pulses is not of type int. :raises ValueError if n_pulses is negative or exceeds the 32bit limit.

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

disable_pulses(**addresses*: Optional[Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel]]) → None

Disable previously configured pulse channels. :param addresses: tuple of FIO addresses. All channels disabled if no argument is passed.

enable_clock(*clock_enabled*: bool) → None

Enable/disable LabJack clock to configure or send pulses. :param clock_enabled: True -> enable, False -> disable. :raises TypeError: if clock_enabled is not of type bool

get_ain(**channels*: int) → Union[float, Sequence[float]]

Read currently measured value (voltage, resistance, ...) from one or more of analog inputs.

Parameters **channels** – AIN number or numbers (0..254)

Returns the read value (voltage, resistance, ...) as float or tuple of them in case multiple channels given

get_cal_current_source(*name*: Union[str, CalMicroAmpere]) → float

This function will return the calibration of the chosen current source, this ist not a measurement!

The value was stored during fabrication.

Parameters **name** – ‘200uA’ or ‘10uA’ current source

Returns calibration of the chosen current source in ampere

get_clock() → Dict[str, Union[int, float]]

Return clock settings read from LabJack.

get_digital_input(*address*: Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel]) → hvl_ccb.dev.labjack.LabJack.DIOStatus

Get the value of a digital input.

allowed names for T7 (Pro): FIO0 - FIO7, EIO0 - EIO 7, CIO0- CIO3, MIO0 - MIO2 :param address: name of the output -> ‘FIO0’ :return: HIGH when address DIO is high, and LOW when address DIO is low

get_product_id() → int

This function returns the product ID reported by the connected device.

Attention: returns 7 for both T7 and T7-Pro devices!

Returns integer product ID of the device

get_product_name(*force_query_id=False*) → str

This function will return the product name based on product ID reported by the device.

Attention: returns “T7” for both T7 and T7-Pro devices!

Parameters `force_query_id` – boolean flag to force `get_product_id` query to device instead of using cached device type from previous queries.

Returns device name string, compatible with `LabJack.DeviceType`

get_product_type(`force_query_id: bool = False`) → `hvl_ccb._dev.labjack.DeviceType`

This function will return the device type based on reported device type and in case of unambiguity based on configuration of device's communication protocol (e.g. for "T7" and "T7_PRO" devices), or, if not available first matching.

Parameters `force_query_id` – boolean flag to force `get_product_id` query to device instead of using cached device type from previous queries.

Returns `DeviceType` instance

Raises `LabJackIdentifierDIOError` – when read Product ID is unknown

get_sbus_rh(`number: int`) → float

Read the relative humidity value from a serial SBUS sensor.

Parameters `number` – port number (0..22)

Returns relative humidity in %RH

get_sbus_temp(`number: int`) → float

Read the temperature value from a serial SBUS sensor.

Parameters `number` – port number (0..22)

Returns temperature in Kelvin

get_serial_number() → int

Returns the serial number of the connected LabJack.

Returns Serial number.

read_resistance(`channel: int`) → float

Read resistance from specified channel.

Parameters `channel` – channel with resistor

Returns resistance value with 2 decimal places

read_thermocouple(`pos_channel: int`) → float

Read the temperature of a connected thermocouple.

Parameters `pos_channel` – is the AIN number of the positive pin

Returns temperature in specified unit

send_pulses(*`addresses: Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel]`) → None

Sends pre-configured pulses for specified addresses. :param `addresses`: tuple of FIO addresses :raises `LabJackError` if an address has not been configured.

set_ain_differential(`pos_channel: int, differential: bool`) → None

Sets an analog input to differential mode or not. T7-specific: For base differential channels, positive must be even channel from 0-12 and negative must be positive+1. For extended channels 16-127, see Mux80 datasheet.

Parameters

- `pos_channel` – is the AIN number (0..12)
- `differential` – True or False

Raises `LabJackError` – if parameters are unsupported

set_ain_range(*channel: int, vrange: Union[Real, AInRange]*) → None

Set the range of an analog input port.

Parameters

- **channel** – is the AIN number (0..254)
- **vrange** – is the voltage range to be set

set_ain_resistance(*channel: int, vrange: Union[Real, AInRange], resolution: int*) → None

Set the specified channel to resistance mode. It utilized the 200uA current source of the LabJack.

Parameters

- **channel** – channel that should measure the resistance
- **vrange** – voltage range of the channel
- **resolution** – resolution index of the channel T4: 0-5, T7: 0-8, T7-Pro 0-12

set_ain_resolution(*channel: int, resolution: int*) → None

Set the resolution index of an analog input port.

Parameters

- **channel** – is the AIN number (0..254)
- **resolution** – is the resolution index within 0...`get_product_type().ain_max_resolution` range; 0 will set the resolution index to default value.

set_ain_thermocouple(*pos_channel: int, thermocouple: Union[None, str, ThermocoupleType], cjc_address: int = 60050, cjc_type: Union[str, CjcType] = CjcType.internal, vrange: Union[Real, AInRange] = AInRange.ONE_HUNDREDTH, resolution: int = 10, unit: Union[str, TemperatureUnit] = TemperatureUnit.K*) → None

Set the analog input channel to thermocouple mode.

Parameters

- **pos_channel** – is the analog input channel of the positive part of the differential pair
- **thermocouple** – None to disable thermocouple mode, or string specifying the thermocouple type
- **cjc_address** – modbus register address to read the CJC temperature
- **cjc_type** – determines cjc slope and offset, ‘internal’ or ‘lm34’
- **vrange** – measurement voltage range
- **resolution** – resolution index (T7-Pro: 0-12)
- **unit** – is the temperature unit to be returned (‘K’, ‘C’ or ‘F’)

Raises `LabJackError` – if parameters are unsupported

set_analog_output(*channel: int, value: Union[int, float]*) → None

Set the voltage of a analog output port

Parameters

- **channel** – DAC channel number 1/0
- **value** – The output voltage value 0-5 Volts int/float

set_clock(*clock_frequency*: Union[Number, ClockFrequency] = 10000000, *clock_period*: Number = 1) → None

Configure LabJack clock for pulse out feature. :param *clock_frequency*: clock frequency in Hz; default 10 MHz for base 10. :raises ValueError: if *clock_frequency* is not allowed (see ClockFrequency). :param *clock_period*: clock roll time in seconds; default 1s, 0 for max. :raises ValueError: if *clock_period* exceeds the 32bit tick limit. Clock period determines pulse spacing when using multi-pulse settings. Ensure period exceeds maximum intended pulse end time.

set_digital_output(*address*: str, *state*: Union[int, DIOStatus]) → None

Set the value of a digital output.

Parameters

- **address** – name of the output -> ‘FIO0’
- **state** – state of the output -> *DIOStatus* instance or corresponding *int* value

start() → None

Start the Device.

stop() → None

Stop the Device.

exception LabJackError

Bases: Exception


Errors of the LabJack device.

exception LabJackIdentifierDIOError

Bases: Exception

Error indicating a wrong DIO identifier

hvl_ccb.dev.lauda

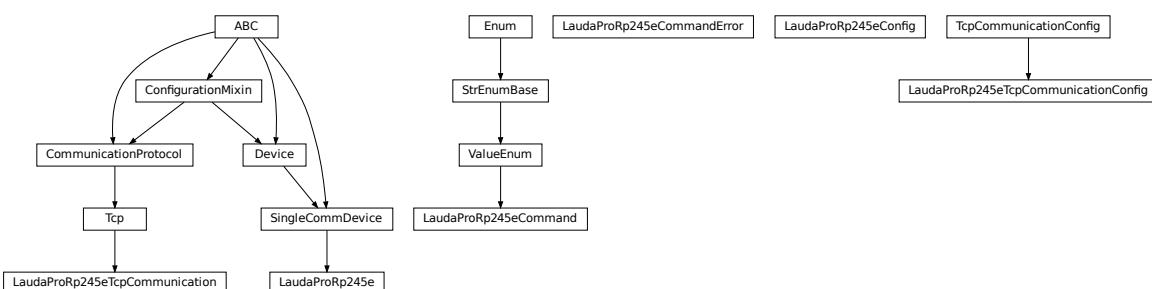

Device class for controlling a Lauda PRO RP245E, circulation chiller over TCP.

class LaudaProRp245e(*com*, *dev_config*=None)

Bases: *hvl_ccb.dev.base.SingleCommDevice*

Lauda RP245E circulation chiller class.

static config_cls() → Type[*hvl_ccb.dev.lauda.LaudaProRp245eConfig*]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

continue_ramp() → str

Continue current ramp program.

Returns reply of the device to the last call of “query”

static default_com_cls() → Type[*hvl_ccb.dev.lauda.LaudaProRp245eTcpCommunication*]

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_bath_temp() → float

:return : float value of measured lauda bath temp in °C

get_device_type() → str

:return : Connected Lauda device type (for connection/com test)

pause() → str

Stop temperature control and pump.

Returns reply of the device to the last call of “query”

pause_ramp() → str

Pause current ramp program.

Returns reply of the device to the last call of “query”

reset_ramp() → str

Delete all segments from current ramp program.

Returns reply of the device to the last call of “query”

run() → str

Start temperature control & pump.

Returns reply of the device to the last call of “query”

set_control_mode(mod: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = ExtControlModeEnum.INTERNAL) → str

Define control mode. 0 = INTERNAL (control bath temp), 1 = EXPT100 (pt100 attached to chiller), 2 = ANALOG, 3 = SERIAL, 4 = USB, 5 = ETH (to be used when passing the ext. temp. via ethernet) (temperature then needs to be passed every second, when not using options 3, 4, or 5)

Parameters **mod** – temp control mode (control internal temp or external temp).

Returns reply of the device to the last call of “query” (“OK”, if command was recognized”)

set_external_temp(external_temp: float = 20.0) → str

Pass value of external controlled temperature. Should be done every second, when control of external temperature is active. Has to be done right before control of external temperature is activated.

Parameters **external_temp** – current value of external temperature to be controlled.

Returns reply of the device to the last call of “query”

set_pump_level(pump_level: int = 6) → str

Set pump level Raises ValueError, if pump level is invalid.

Parameters **pump_level** – pump level.

Returns reply of the device to the last call of “query”

set_ramp_iterations(num: int = 1) → str

Define number of ramp program cycles.

Parameters **num** – number of program cycles to be performed.

Returns reply of the device to the last call of “query”

set_ramp_program(*program: int = 1*) → str

Define ramp program for following ramp commands. Raises ValueError if maximum number of ramp programs (5) is exceeded.

Parameters **program** – Number of ramp program to be activated for following commands.

Returns reply of the device to the last call of “query”

set_ramp_segment(*temp: float = 20.0, dur: int = 0, tol: float = 0.0, pump: int = 6*) → str

Define segment of current ramp program - will be attached to current program. Raises ValueError, if pump level is invalid.

Parameters

- **temp** – target temperature of current ramp segment
- **dur** – duration in minutes, in which target temperature should be reached
- **tol** – tolerance at which target temperature should be reached (for 0.00, next segment is started after dur has passed).
- **pump** – pump level to be used for this program segment.

Returns reply of the device to the last call of “query”

set_temp_set_point(*temp_set_point: float = 20.0*) → str

Define temperature set point

Parameters **temp_set_point** – temperature set point.

Returns reply of the device to the last call of “query”

start() → None

Start this device.

start_ramp() → str

Start current ramp program.

Returns reply of the device to the last call of “query”

stop() → None

Stop this device. Disables access and closes the communication protocol.

stop_ramp() → str

Stop current ramp program.

Returns reply of the device to the last call of “query”

validate_pump_level(*level: int*)

Validates pump level. Raises ValueError, if pump level is incorrect. :param level: pump level, integer

class LaudaProRp245eCommand(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: [hvl_ccb.utils.enum.ValueEnum](#)

Commands for Lauda PRO RP245E Chiller Command strings most often need to be complimented with a parameter (attached as a string) before being sent to the device. Commands implemented as defined in “Lauda Betriebsanleitung fuer PRO Badthermostate und Umwaelzthermostate” pages 42 - 49

BATH_TEMP = 'IN_PV_00'

Request internal bath temperature

COM_TIME_OUT = 'OUT_SP_08_'

Define communication time out

```

CONT_MODE = 'OUT_MODE_01_'
    Set control mode 1=internal, 2=ext. analog, 3=ext. serial, 4=USB, 5=ethernet

DEVICE_TYPE = 'TYPE'
    Request device type

EXTERNAL_TEMP = 'OUT_PV_05_'
    Pass on external controlled temperature

LOWER_TEMP = 'OUT_SP_05_'
    Define lower temp limit

OPERATION_MODE = 'OUT_SP_02_'
    Define operation mode

PUMP_LEVEL = 'OUT_SP_01_'
    Define pump level 1-8

RAMP_CONTINUE = 'RMP_CONT'
    Continue a paused ramp program

RAMP_DELETE = 'RMP_RESET'
    Reset a selected ramp program

RAMP_ITERATIONS = 'RMP_OUT_02_'
    Define how often a ramp program should be iterated

RAMP_PAUSE = 'RMP_PAUSE'
    Pause a selected ramp program

RAMP_SELECT = 'RMP_SELECT_'
    Select a ramp program (target for all further ramp commands)

RAMP_SET = 'RMP_OUT_00_'
    Define parameters of a selected ramp program

RAMP_START = 'RMP_START'
    Start a selected ramp program

RAMP_STOP = 'RMP_STOP'
    Stop a running ramp program

START = 'START'
    Start temp control (pump and heating/cooling)

STOP = 'STOP'
    Stop temp control (pump and heating/cooling)

TEMP_SET_POINT = 'OUT_SP_00_'
    Define temperature set point

UPPER_TEMP = 'OUT_SP_04_'
    Define upper temp limit

build_str(param: str = "", terminator: str = '\r\n')
    Build a command string for sending to the device

```

Parameters

- **param** – Command's parameter given as string
- **terminator** – Command's terminator

Returns Command's string with a parameter and terminator

```
exception LaudaProRp245eCommandError
```

Bases: Exception

Exception raised when an error is returned upon a command.

```
class LaudaProRp245eConfig(temp_set_point_init: Union[int, float] = 20.0, pump_init: int = 6, upper_temp: Union[int, float] = 80.0, lower_temp: Union[int, float] = - 55.0, com_time_out: Union[int, float] = 0, max_pump_level: int = 8, max_pr_number: int = 5, operation_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.OperationModeEnum] = OperationModeEnum.AUTO, control_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = ExtControlModeEnum.INTERNAL)
```

Bases: object

Configuration for the Lauda RP245E circulation chiller.

```
class ExtControlModeEnum(value)
```

Bases: enum.IntEnum

Source for definition of external, controlled temperature (option 2, 3 and 4 are not available with current configuration of the Lauda RP245E, add-on hardware would required)

```
ANALOG = 2
```

```
ETH = 5
```

```
EXPT100 = 1
```

```
INTERNAL = 0
```

```
SERIAL = 3
```

```
USB = 4
```

```
class OperationModeEnum(value)
```

Bases: enum.IntEnum

Operation Mode (Cooling OFF/Cooling On/AUTO - set to AUTO)

```
AUTO = 2
```

Automatically select heating/cooling

```
COOLOFF = 0
```

```
COOLON = 1
```

```
clean_values() → None
```

```
com_time_out: Union[int, float] = 0
```

Communication time out (0 = OFF)

```
control_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = 0
```

```
force_value(fieldname, value)
```

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
is_configdataclass = True

classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

lower_temp: Union[int, float] = -55.0
    Lower temperature limit (safe for Galden HT135 cooling liquid)

max_pr_number: int = 5
    Maximum number of ramp programs that can be stored in the memory of the chiller

max_pump_level: int = 8
    Highest pump level of the chiller

operation_mode: Union[int,
hvl_ccb.dev.lauda.LaudaProRp245eConfig.OperationModeEnum] = 2

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

pump_init: int = 6
    Default pump Level

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

temp_set_point_init: Union[int, float] = 20.0
    Default temperature set point

upper_temp: Union[int, float] = 80.0
    Upper temperature limit (safe for Galden HT135 cooling liquid)

class LaudaProRp245eTcpCommunication(configuration)
    Bases: hvl\_ccb.comm.tcp.Tcp

    Implements the Communication Protocol for Lauda PRO RP245E TCP connection.

close() → None
    Close the Lauda PRO RP245E TCP connection.

static config_cls() → Type[hvl\_ccb.dev.lauda.LaudaProRp245eTcpCommunicationConfig]
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

open() → None
    Open the Lauda PRO RP245E TCP connection.

    Raises LaudaProRp245eCommandError – if the connection fails.

query_command(command: hvl_ccb.dev.lauda.LaudaProRp245eCommand, param: str = "") → str
    Send and receive function. E.g. to be used when setting/changing device setting. :param command: first part of command string, defined in LaudaProRp245eCommand :param param: second part of command string, parameter (by default '') :return: None
```

read() → str

Receive value function. :return: reply from device as a string, the terminator, as well as the ‘OK’ stripped from the reply to make it directly useful as a value (e.g. in case the internal bath temperature is requested)

write_command(command: hvl_ccb.dev.lauda.LaudaProRp245eCommand, param: str = "") → None

Send command function. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None

class LaudaProRp245eTcpCommunicationConfig(host: str, port: int = 54321, bufsize: int = 1024, wait_sec_pre_read_or_write: Union[int, float] = 0.005, terminator: str = '\r\n')

Bases: *hvl_ccb.comm.tcp.TcpCommunicationConfig*

Configuration dataclass for *LaudaProRp245eTcpCommunication*.

clean_values() → None**force_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns a list of strings containing all required keys.

terminator: str = '\r\n'

The terminator character

wait_sec_pre_read_or_write: Union[int, float] = 0.005

Delay time between commands in seconds

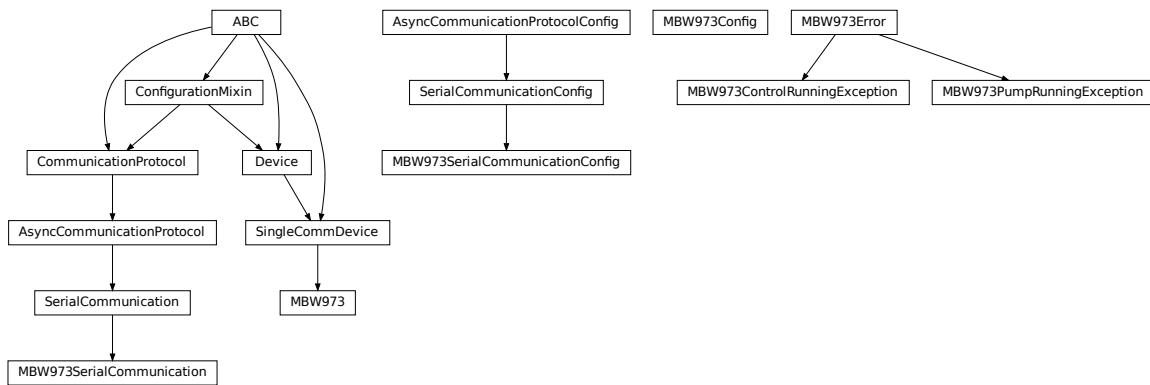hvl_ccb.dev.mbw973

Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

class MBW973(com, dev_config=None)

Bases: `hvl_ccb.dev.base.SingleCommDevice`

MBW 973 dew point mirror device class.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

is_done() → bool

Poll status of the dew point mirror and return True, if all measurements are done.

Returns True, if all measurements are done; False otherwise.

Raises `SerialCommunicationIOError` – when communication port is not opened

read(cast_type: Type = <class 'str'>)

Read value from `self.com` and cast to `cast_type`. Raises `ValueError` if read text (`str`) is not convertible to `cast_type`, e.g. to `float` or to `int`.

Returns Read value of `cast_type` type.

read_float() → float

Convenience wrapper for `self.read()`, with typing hint for return value.

Returns Read `float` value.

read_int() → int

Convenience wrapper for `self.read()`, with typing hint for return value.

Returns Read `int` value.

read_measurements() → Dict[str, float]

Read out measurement values and return them as a dictionary.

Returns Dictionary with values.

Raises *SerialCommunicationIOError* – when communication port is not opened

set_measuring_options(humidity: bool = True, sf6_purity: bool = False) → None

Send measuring options to the dew point mirror.

Parameters

- **humidity** – Perform humidity test or not?
- **sf6_purity** – Perform SF6 purity test or not?

Raises *SerialCommunicationIOError* – when communication port is not opened

start() → None

Start this device. Opens the communication protocol and retrieves the set measurement options from the device.

Raises *SerialCommunicationIOError* – when communication port cannot be opened.

start_control() → None

Start dew point control to acquire a new value set.

Raises *SerialCommunicationIOError* – when communication port is not opened

stop() → None

Stop the device. Closes also the communication protocol.

write(value) → None

Send *value* to *self.com*.

Parameters **value** – Value to send, converted to *str*.

Raises *SerialCommunicationIOError* – when communication port is not opened

class MBW973Config(polling_interval: Union[int, float] = 2)

Bases: object

Device configuration dataclass for MBW973.

clean_values()

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

polling_interval: Union[int, float] = 2
 Polling period for *is_done* status queries [in seconds].

classmethod required_keys() → Sequence[str]
 Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception MBW973ControlRunningException
 Bases: *hvl_ccb.dev.mbw973.MBW973Error*
 Error indicating there is still a measurement running, and a new one cannot be started.

exception MBW973Error
 Bases: *Exception*
 General error with the MBW973 dew point mirror device.

exception MBW973PumpRunningException
 Bases: *hvl_ccb.dev.mbw973.MBW973Error*
 Error indicating the pump of the dew point mirror is still recovering gas, unable to start a new measurement.

class MBW973SerialCommunication(configuration)
 Bases: *hvl_ccb.comm.serial.SerialCommunication*
 Specific communication protocol implementation for the MBW973 dew point mirror. Already predefines device-specific protocol parameters in config.

static config_cls()
 Return the default configdataclass class.

Returns a reference to the default configdataclass class

class MBW973SerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)
 Bases: *hvl_ccb.comm.serial.SerialCommunicationConfig*

baudrate: int = 9600
 Baudrate for MBW973 is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
 One byte is eight bits long

force_value(fieldname, value)
 Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, [hvl_ccb.comm.serial.SerialCommunicationParity](#)] = 'N'

MBW973 does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, [hvl_ccb.comm.serial.SerialCommunicationStopbits](#)] = 1

MBW973 does use one stop bit


terminator: bytes = b'\r'

The terminator is only CR

timeout: Union[int, float] = 3

use 3 seconds timeout as default

[hvl_ccb.dev.newport](#)

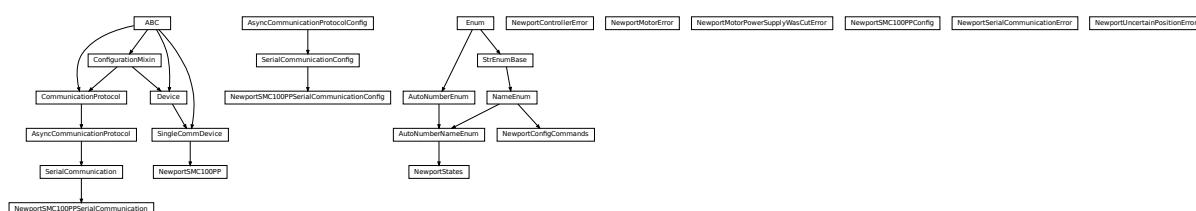

Device class for Newport SMC100PP stepper motor controller with serial communication.

The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/f/sm100-single-axis-dc-or-stepper-motion-controller>

class NewportConfigCommands(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [hvl_ccb.utils.enum.NameEnum](#)

Commands predefined by the communication protocol of the SMC100PP

AC = 'acceleration'

```

BA = 'backlash_compensation'
BH = 'hysteresis_compensation'
FRM = 'micro_step_per_full_step_factor'
FRS = 'motion_distance_per_full_step'
HT = 'home_search_type'
JR = 'jerk_time'
OH = 'home_search_velocity'
OT = 'home_search_timeout'
QIL = 'peak_output_current_limit'
SA = 'rs485_address'
SL = 'negative_software_limit'
SR = 'positive_software_limit'
VA = 'velocity'
VB = 'base_velocity'
ZX = 'stage_configuration'

exception NewportControllerError
    Bases: Exception
    Error with the Newport controller.

exception NewportMotorError
    Bases: Exception
    Error with the Newport motor.

exception NewportMotorPowerSupplyWasCutError
    Bases: Exception
    Error with the Newport motor after the power supply was cut and then restored, without interrupting the communication with the controller.

class NewportSMC100PP(com, dev_config=None)
    Bases: hvl\_ccb.dev.base.SingleCommDevice
    Device class of the Newport motor controller SMC100PP

class MotorErrors(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: aenum.Enum
    Possible motor errors reported by the motor during get_state().
    DC_VOLTAGE_TOO_LOW = 3
    FOLLOWING_ERROR = 6
    HOMING_TIMEOUT = 5
    NED_END_OF_TURN = 11
    OUTPUT_POWER_EXCEEDED = 2
    PEAK_CURRENT_LIMIT = 9
    POS_END_OF_TURN = 10

```

```
RMS_CURRENT_LIMIT = 8
SHORT_CIRCUIT = 7
WRONG_ESP_STAGE = 4

class StateMessages(value=<no_arg>, names=None, module=None, type=None, start=1,
                    boundary=None)
Bases: aenum.Enum

Possible messages returned by the controller on get_state() query.

CONFIG = '14'
DISABLE_FROM_JOGGING = '3E'
DISABLE_FROM_MOVING = '3D'
DISABLE_FROM_READY = '3C'
HOMING_FROM_RS232 = '1E'
HOMING_FROM_SMC = '1F'
JOGGING_FROM_DISABLE = '47'
JOGGING_FROM_READY = '46'
MOVING = '28'
NO_REF_ESP_STAGE_ERROR = '10'
NO_REF_FROM_CONFIG = '0C'
NO_REF_FROM_DISABLED = '0D'
NO_REF_FROM_HOMING = '0B'
NO_REF_FROM_JOGGING = '11'
NO_REF_FROM_MOVING = '0F'
NO_REF_FROM_READY = '0E'
NO_REF_FROM_RESET = '0A'
READY_FROM_DISABLE = '34'
READY_FROM_HOMING = '32'
READY_FROM_JOGGING = '35'
READY_FROM_MOVING = '33'

States
alias of hvl\_ccb.dev.newport.NewportStates

static config_cls()
Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device
```

exit_configuration(*add*: *Optional[int]* = *None*) → *None*

Exit the CONFIGURATION state and go back to the NOT REFERENCED state. All configuration parameters are saved to the device's memory.

Parameters **add** – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_acceleration(*add*: *Optional[int]* = *None*) → *Union[int, float]*

Leave the configuration state. The configuration parameters are saved to the device's memory.

Parameters **add** – controller address (1 to 31)

Returns acceleration (preset units/s^2), value between 1e-6 and 1e12

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_controller_information(*add*: *Optional[int]* = *None*) → *str*

Get information on the controller name and driver version

Parameters **add** – controller address (1 to 31)

Returns controller information

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_motor_configuration(*add*: *Optional[int]* = *None*) → *Dict[str, float]*

Query the motor configuration and returns it in a dictionary.

Parameters **add** – controller address (1 to 31)

Returns dictionary containing the motor's configuration

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_move_duration(*dist*: *Union[int, float]*, *add*: *Optional[int]* = *None*) → *float*

Estimate the time necessary to move the motor of the specified distance.

Parameters

- **dist** – distance to travel
- **add** – controller address (1 to 31), defaults to self.address

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_negative_software_limit(*add: Optional[int] = None*) → Union[int, float]

Get the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

Parameters *add* – controller address (1 to 31)

Returns negative software limit (preset units), value between -1e12 and 0

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_position(*add: Optional[int] = None*) → float

Returns the value of the current position.

Parameters *add* – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportUncertainPositionError* – if the position is ambiguous

get_positive_software_limit(*add: Optional[int] = None*) → Union[int, float]

Get the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

Parameters *add* – controller address (1 to 31)

Returns positive software limit (preset units), value between 0 and 1e12

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_state(*add: int = None*) → *StateMessages*

Check on the motor errors and the controller state

Parameters *add* – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportMotorError* – if the motor reports an error

Returns state message from the device (member of StateMessages)

`go_home(add: Optional[int] = None) → None`

Move the motor to its home position.

Parameters `add` – controller address (1 to 31), defaults to self.address

Raises

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

`go_to_configuration(add: Optional[int] = None) → None`

This method is executed during start(). It can also be executed after a reset(). The controller is put in CONFIG state, where configuration parameters can be changed.

Parameters `add` – controller address (1 to 31)

Raises

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

`initialize(add: Optional[int] = None) → None`

Puts the controller from the NOT_REF state to the READY state. Sends the motor to its “home” position.

Parameters `add` – controller address (1 to 31)

Raises

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

`move_to_absolute_position(pos: Union[int, float], add: Optional[int] = None) → None`

Move the motor to the specified position.

Parameters

- `pos` – target absolute position (affected by the configured offset)
- `add` – controller address (1 to 31), defaults to self.address

Raises

- `SerialCommunicationIOError` – if the com is closed
- `NewportSerialCommunicationError` – if an unexpected answer is obtained
- `NewportControllerError` – if the controller reports an error

`move_to_relative_position(pos: Union[int, float], add: Optional[int] = None) → None`

Move the motor of the specified distance.

Parameters

- `pos` – distance to travel (the sign gives the direction)
- `add` – controller address (1 to 31), defaults to self.address

Raises

- `SerialCommunicationIOError` – if the com is closed

- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

reset(*add: Optional[int] = None*) → None

Resets the controller, equivalent to a power-up. This puts the controller back to NOT REFERENCED state, which is necessary for configuring the controller.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_acceleration(*acc: Union[int, float], add: Optional[int] = None*) → None

Leave the configuration state. The configuration parameters are saved to the device's memory.

Parameters

- **acc** – acceleration (preset units/s^2), value between 1e-6 and 1e12
- **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_motor_configuration(*add: Optional[int] = None, config: Optional[dict] = None*) → None

Set the motor configuration. The motor must be in CONFIG state.

Parameters

- **add** – controller address (1 to 31)
- **config** – dictionary containing the motor's configuration

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_negative_software_limit(*lim: Union[int, float], add: Optional[int] = None*) → None

Set the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

Parameters

- **lim** – negative software limit (preset units), value between -1e12 and 0
- **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_positive_software_limit(*lim: Union[int, float]*, *add: Optional[int] = None*) → None

Set the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

Parameters

- **lim** – positive software limit (preset units), value between 0 and 1e12
- **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

start()

Opens the communication protocol and applies the config.

Raises **SerialCommunicationIOError** – when communication port cannot be opened

stop() → None

Stop the device. Close the communication protocol.

stop_motion(*add: Optional[int] = None*) → None

Stop a move in progress by decelerating the positioner immediately with the configured acceleration until it stops. If a controller address is provided, stops a move in progress on this controller, else stops the moves on all controllers.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

wait_until_motor_initialized(*add: Optional[int] = None*) → None

Wait until the motor leaves the HOMING state (at which point it should have arrived to the home position).

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

```
class NewportSMC100PPConfig(address: int = 1, user_position_offset: Union[int, float] = 23.987,
    screw_scaling: Union[int, float] = 1, exit_configuration_wait_sec: Union[int,
    float] = 5, move_wait_sec: Union[int, float] = 1, acceleration: Union[int, float]
    = 10, backlash_compensation: Union[int, float] = 0, hysteresis_compensation:
    Union[int, float] = 0.015, micro_step_per_full_step_factor: int = 100,
    motion_distance_per_full_step: Union[int, float] = 0.01, home_search_type:
    Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] =
    HomeSearch.HomeSwitch, jerk_time: Union[int, float] = 0.04,
    home_search_velocity: Union[int, float] = 4, home_search_timeout: Union[int,
    float] = 27.5, home_search_polling_interval: Union[int, float] = 1,
    peak_output_current_limit: Union[int, float] = 0.4, rs485_address: int = 2,
    negative_software_limit: Union[int, float] = - 23.5, positive_software_limit:
    Union[int, float] = 25, velocity: Union[int, float] = 4, base_velocity: Union[int,
    float] = 0, stage_configuration: Union[int,
    hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageConfig] =
    EspStageConfig.EnableEspStageCheck)
```

Bases: `object`

Configuration dataclass for the Newport motor controller SMC100PP.

```
class EspStageConfig(value=<no_arg>, names=None, module=None, type=None, start=1,
    boundary=None)
```

Bases: `aenum.IntEnum`

Different configurations to check or not the motor configuration upon power-up.

`DisableEspStageCheck` = 1

`EnableEspStageCheck` = 3

`UpdateEspStageInfo` = 2

```
class HomeSearch(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `aenum.IntEnum`

Different methods for the motor to search its home position during initialization.

`CurrentPosition` = 1

`EndOfRunSwitch` = 4

`EndOfRunSwitch_and_Index` = 3

`HomeSwitch` = 2

`HomeSwitch_and_Index` = 0

`acceleration: Union[int, float]` = 10

`address: int` = 1

`backlash_compensation: Union[int, float]` = 0

`base_velocity: Union[int, float]` = 0

`clean_values()`

`exit_configuration_wait_sec: Union[int, float]` = 5

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

home_search_polling_interval: Union[int, float] = 1

home_search_timeout: Union[int, float] = 27.5

home_search_type: Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] = 2

home_search_velocity: Union[int, float] = 4

hysteresis_compensation: Union[int, float] = 0.015

is_configdataclass = True

jerk_time: Union[int, float] = 0.04

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

micro_step_per_full_step_factor: int = 100

motion_distance_per_full_step: Union[int, float] = 0.01

property motor_config: Dict[str, float]

Gather the configuration parameters of the motor into a dictionary.

Returns dict containing the configuration parameters of the motor

move_wait_sec: Union[int, float] = 1

negative_software_limit: Union[int, float] = -23.5

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

peak_output_current_limit: Union[int, float] = 0.4

positive_software_limit: Union[int, float] = 25

post_force_value(fieldname, value)

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

rs485_address: int = 2

screw_scaling: Union[int, float] = 1

stage_configuration: Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageConfig] = 3

user_position_offset: Union[int, float] = 23.987

velocity: Union[int, float] = 4

```
class NewportSMC100PPSerialCommunication(configuration)
```

Bases: [hvl_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for NewportSMC100 controller. Already predefines device-specific protocol parameters in config.

```
class ControllerErrors(value=<no_arg>, names=None, module=None, type=None, start=1,  
boundary=None)
```

Bases: [aenum.Enum](#)

Possible controller errors with values as returned by the device in response to sent commands.

ADDR_INCORRECT = 'B'

CMD_EXEC_ERROR = 'V'

CMD_NOT_ALLOWED = 'D'

CMD_NOT_ALLOWED_CC = 'X'

CMD_NOT_ALLOWED_CONFIGURATION = 'I'

CMD_NOT_ALLOWED_DISABLE = 'J'

CMD_NOT_ALLOWED_HOMING = 'L'

CMD_NOT_ALLOWED_MOVING = 'M'

CMD_NOT_ALLOWED_NOT_REFERENCED = 'H'

CMD_NOT_ALLOWED_PP = 'W'

CMD_NOT_ALLOWED_READY = 'K'

CODE_OR_ADDR_INVALID = 'A'

COM_TIMEOUT = 'S'

DISPLACEMENT_OUT_OF_LIMIT = 'G'

EEPROM_ACCESS_ERROR = 'U'

ESP_STAGE_NAME_INVALID = 'F'

HOME_STARTED = 'E'

NO_ERROR = '@'

PARAM_MISSING_OR_INVALID = 'C'

POSITION_OUT_OF_LIMIT = 'N'

check_for_error(add: int) → None

Ask the Newport controller for the last error it recorded.

This method is called after every command or query.

Parameters add – controller address (1 to 31)

Raises

- [SerialCommunicationIOError](#) – if the com is closed
- [NewportSerialCommunicationError](#) – if an unexpected answer is obtained
- [NewportControllerError](#) – if the controller reports an error

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(*add*: int, *cmd*: str, *param*: Optional[Union[int, float, str]] = None) → str

Send a query to the controller, read the answer, and check for errors. The prefix add+cmd is removed from the answer.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

Returns the answer from the device without the prefix

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

query_multiple(*add*: int, *cmd*: str, *prefixes*: List[str]) → List[str]

Send a query to the controller, read the answers, and check for errors. The prefixes are removed from the answers.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **prefixes** – prefixes of each line expected in the answer

Returns list of answers from the device without prefix

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

read_text() → str

Read one line of text from the serial port, and check for presence of a null char which indicates that the motor power supply was cut and then restored. The input buffer may hold additional data afterwards, since only one line is read.

This method uses *self.access_lock* to ensure thread-safety.

Returns String read from the serial port; '' if there was nothing to read.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **NewportMotorPowerSupplyWasCutError** – if a null char is read

send_command(*add*: int, *cmd*: str, *param*: Optional[Union[int, float, str]] = None) → None

Send a command to the controller, and check for errors.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent

- **param** – optional parameter (int/float/str) appended to the command

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

send_stop(*add*: int) → None

Send the general stop ST command to the controller, and check for errors.

Parameters *add* – the controller address (1 to 31)**Returns** ControllerErrors reported by Newport Controller**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained

```
class NewportSMC100PPSSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'ascii',
                                                encoding_error_handling: str = 'replace',
                                                wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                                default_n_attempts_read_text_nonempty: int = 10,
                                                port: Union[str, NoneType] = None, baudrate: int =
                                                57600, parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 10)
```

Bases: *hvl_ccb.comm.serial.SerialCommunicationConfig***baudrate: int = 57600**

Baudrate for NewportSMC100 controller is 57600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

NewportSMC100 controller uses 8 bits for one data byte

encoding: str = 'ascii'

use ASCII as de-/encoding, cf. the manual

encoding_error_handling: str = 'replace'

replace bytes with instead of raising utf-8 exception when decoding fails

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

```
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'
    NewportSMC100 controller does not use parity

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
    NewportSMC100 controller uses one stop bit

terminator: bytes = b'\r\n'
    The terminator is CR/LF

timeout: Union[int, float] = 10
    use 10 seconds timeout as default

exception NewportSerialCommunicationError
    Bases: Exception

    Communication error with the Newport controller.

class NewportStates(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl_ccb.utils.enum.AutoNumberNameEnum


    States of the Newport controller. Certain commands are allowed only in certain states.

    CONFIG = 3
    DISABLE = 6
    HOMING = 2
    JOGGING = 7
    MOVING = 5
    NO_REF = 1
    READY = 4

exception NewportUncertainPositionError
    Bases: Exception

    Error with the position of the Newport motor.
```

`hvl_ccb.dev.pfeiffer_tpg`

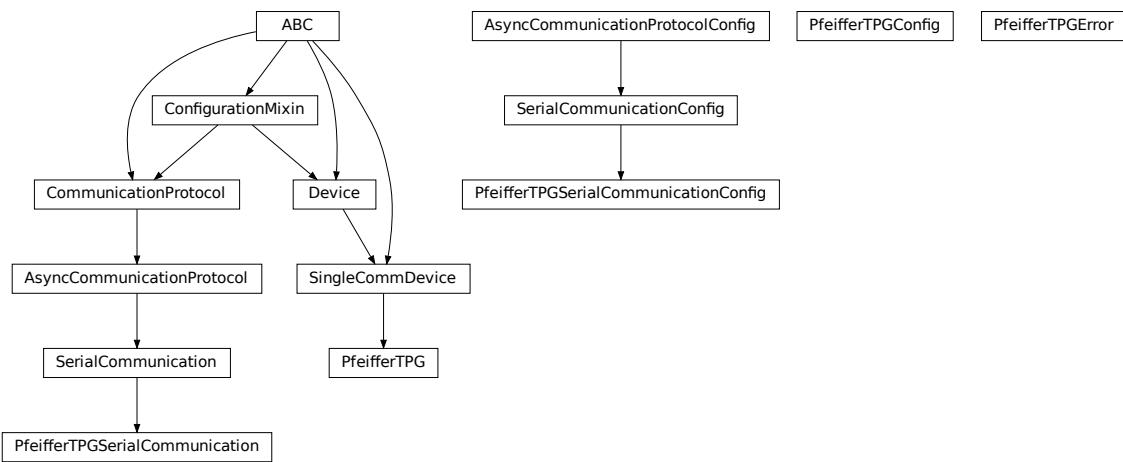

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/> measurement/activeline/controllers/

class PfeifferTPG(*com, dev_config=None*)
Bases: *hvl_ccb.dev.base.SingleCommDevice*

Pfeiffer TPG control unit device class

class PressureUnits(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)
Bases: *hvl_ccb.utils.enum.NameEnum*

Enum of available pressure units for the digital display. “0” corresponds either to bar or to mbar depending on the TPG model. In case of doubt, the unit is visible on the digital display.

Micron = 3

Pascal = 2

Torr = 1

Volt = 5

bar = 0

hPascal = 4

mbar = 0

class SensorStatus(*value*)

Bases: *enum.IntEnum*

An enumeration.

Identification_error = 6

No_sensor = 5

```

Ok = 0
OVERRANGE = 2
Sensor_error = 3
Sensor_off = 4
Underrange = 1

class SensorTypes(value)
    Bases: enum.Enum

    An enumeration.

    CMR = 4
    IKR = 2
    IKR11 = 2
    IKR9 = 2
    IMR = 5
    None = 7
    PBR = 6
    PKR = 3
    TPR = 1
    noSENSOR = 7
    noSen = 7

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

get_full_scale_mbar() → List[Union[int, float]]
    Get the full scale range of the attached sensors

    Returns full scale range values in mbar, like [0.01, 1, 0.1, 1000, 50000, 10]

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • PfeifferTPGError – if command fails

get_full_scale_unitless() → List[int]
    Get the full scale range of the attached sensors. See lookup table between command and corresponding
    pressure in the device user manual.

    Returns list of full scale range values, like [0, 1, 3, 3, 2, 0]

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • PfeifferTPGError – if command fails

```

identify_sensors() → None

Send identification request TID to sensors on all channels.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

measure(channel: int) → Tuple[str, float]

Get the status and measurement of one sensor

Parameters **channel** – int channel on which the sensor is connected, with $1 \leq \text{channel} \leq \text{number_of_sensors}$

Returns measured value as float if measurement successful, sensor status as string if not

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

measure_all() → List[Tuple[str, float]]

Get the status and measurement of all sensors (this command is not available on all models)

Returns list of measured values as float if measurements successful, and or sensor status as strings if not

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

property number_of_sensors

set_display_unit(unit: Union[str, hvl_ccb.dev.pfeiffer_tpg.PfeifferTPG.PressureUnits]) → None

Set the unit in which the measurements are shown on the display.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

set_full_scale_mbar(fsr: List[Union[int, float]]) → None

Set the full scale range of the attached sensors (in unit mbar)

Parameters **fsr** – full scale range values in mbar, for example [0.01, 1000]

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

set_full_scale_unitless(fsr: List[int]) → None

Set the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

Parameters **fsr** – list of full scale range values, like [0, 1, 3, 3, 2, 0]

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

start() → None

Start this device. Opens the communication protocol, and identify the sensors.

Raises `SerialCommunicationIOError` – when communication port cannot be opened

stop() → None

Stop the device. Closes also the communication protocol.

property unit

The pressure unit of readings is always mbar, regardless of the display unit.

class PfeifferTPGConfig(*model*: Union[str, `hvl_ccb.dev.pfeiffer_tpg.PfeifferTPGConfig.Model`] = `Model.TPG25xA`)

Bases: object

Device configuration dataclass for Pfeiffer TPG controllers.

class Model(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: `hvl_ccb.utils.enum.NameEnum`

An enumeration.

`TPG25xA` = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

`TPGx6x` = {0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8, 50000: 9}

is_valid_scale_range_reversed_str(*v*: str) → bool

Check if given string represents a valid reversed scale range of a model.

Parameters *v* – Reversed scale range string.

Returns *True* if valid, *False* otherwise.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

model: Union[str, `hvl_ccb.dev.pfeiffer_tpg.PfeifferTPGConfig.Model`] = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception PfeifferTPGError

Bases: Exception

Error with the Pfeiffer TPG Controller.

class PfeifferTPGSerialCommunication(configuration)

Bases: *hvl_ccb.comm.serial.SerialCommunication*

Specific communication protocol implementation for Pfeiffer TPG controllers. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(cmd: str) → str

Send a query, then read and returns the first line from the com port.

Parameters cmd – query message to send to the device

Returns first line read on the com

Raises

- *SerialCommunicationIOError* – when communication port is not opened

- *PfeifferTPGError* – if the device does not acknowledge the command or if the answer from the device is empty

send_command(cmd: str) → None

Send a command to the device and check for acknowledgement.

Parameters cmd – command to send to the device

Raises

- *SerialCommunicationIOError* – when communication port is not opened

- *PfeifferTPGError* – if the answer from the device differs from the expected acknowledgement character ‘chr(6)’.

```
class PfeifferTPGSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                             Union[str,
                                               hvl_ccb.comm.serial.SerialCommunicationParity] =
                                               <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                               Union[int,
                                                 hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                 <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                               Union[int,
                                                 hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                 <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                               Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for Pfeiffer TPG controllers is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

Pfeiffer TPG controllers do not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1

Pfeiffer TPG controllers use one stop bit

terminator: bytes = b'\r\n'

The terminator is <CR><LF>

```
timeout: Union[int, float] = 3
use 3 seconds timeout as default
```

hvl_ccb.dev.picotech_pt104


NOTE: PicoSDK Python wrappers already on import attempt to load the PicoSDK library; thus, the API docs can only be generated in a system with the latter installed and are by default disabled.

To build the API docs for this submodule locally edit the `docs/hvl_ccb.dev.picotech_pt104.rst` file to remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.picotech_pt104
:parts: 1

.. automodule:: hvl_ccb.dev.picotech_pt104
:members:
:undoc-members:
:show-inheritance:
```

hvl_ccb.dev.rs_rto1024

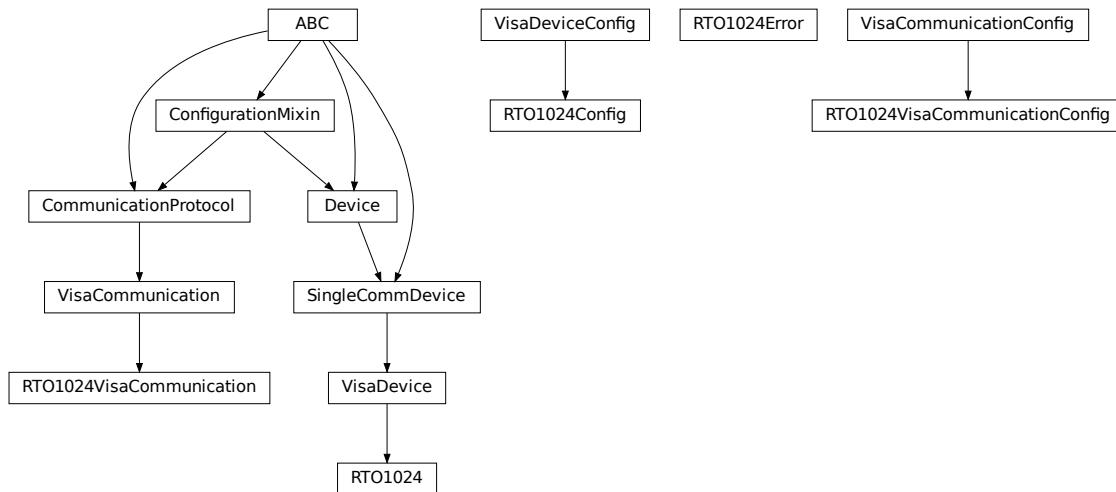

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

```
class RT01024(com: Union[hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunication,
                        hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunicationConfig, dict], dev_config:
                        Union[hvl_ccb.dev.rs_rto1024.RTO1024Config, dict])
Bases: hvl_ccb.dev.visa.VisaDevice

Device class for the Rhode & Schwarz RTO 1024 oscilloscope.
```

```
class TriggerModes(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl\_ccb.utils.enum.AutoNumberNameEnum

    Enumeration for the three available trigger modes.

    AUTO = 1
    FREEERUN = 3
    NORMAL = 2

    classmethod names()
        Returns a list of the available trigger modes. :return: list of strings

activate_measurements(meas_n: int, source: str, measurements: List[str], category: str = 'AMPTime')
    Activate the list of ‘measurements’ of the waveform ‘source’ in the measurement box number ‘meas_n’. The list ‘measurements’ starts with the main measurement and continues with additional measurements of the same ‘category’.

Parameters

- meas_n – measurement number 1..8
- source – measurement source, for example C1W1
- measurements – list of measurements, the first one will be the main measurement.
- category – the category of measurements, by default AMPTime

backup_waveform(filename: str) → None
    Backup a waveform file from the standard directory specified in the device configuration to the standard backup destination specified in the device configuration. The filename has to be specified without .bin or path.

Parameters filename – The waveform filename without extension and path

static config_cls()
    Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
    Return the default communication protocol for this device type, which is VisaCommunication.

Returns the VisaCommunication class

file_copy(source: str, destination: str) → None
    Copy a file from one destination to another on the oscilloscope drive. If the destination file already exists, it is overwritten without notice.

Parameters

- source – absolute path to the source file on the DSO filesystem
- destination – absolute path to the destination file on the DSO filesystem

Raises RT01024Error – if the operation did not complete

get_acquire_length() → float
    Gets the time of one acquisition, that is the time across the 10 divisions of the diagram.



- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]

```

Returns the time for one acquisition. Range: 250e-12 ... 500 [s]

get_channel_offset(*channel: int*) → float

Gets the voltage offset of the indicated channel.

Parameters **channel** – is the channel number (1..4)

Returns channel offset voltage in V (value between -1 and 1)

get_channel_position(*channel: int*) → float

Gets the vertical position of the indicated channel.

Parameters **channel** – is the channel number (1..4)

Returns channel position in div (value between -5 and 5)

get_channel_range(*channel: int*) → float

Queries the channel range in V.

Parameters **channel** – is the input channel (1..4)

Returns channel range in V

get_channel_scale(*channel: int*) → float

Queries the channel scale in V/div.

Parameters **channel** – is the input channel (1..4)

Returns channel scale in V/div

get_channel_state(*channel: int*) → bool

Queries if the channel is active or not.

Parameters **channel** – is the input channel (1..4)

Returns True if active, else False

get_reference_point() → int

Gets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 ... 100 [%]
- Increment: 1 [%]

Returns the reference in %

get_repetitions() → int

Get the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 ... 16777215
- Increment: 10
- *RST = 1

Returns the number of waveforms to acquire

get_timestamps() → List[float]

Gets the timestamps of all recorded frames in the history and returns them as a list of floats.

Returns list of timestamps in [s]

Raises **RT01024Error** – if the timestamps are invalid

list_directory(*path: str*) → List[Tuple[str, str, int]]

List the contents of a given directory on the oscilloscope filesystem.

Parameters **path** – is the path to a folder

Returns a list of filenames in the given folder

load_configuration(*filename: str*) → None

Load current settings from a configuration file. The filename has to be specified without base directory and ‘.dfl’ extension.

Information from the manual *ReCaLl* calls up the instrument settings from an intermediate memory identified by the specified number. The instrument settings can be stored to this memory using the command *\$AV with the associated number. It also activates the instrument settings which are stored in a file and loaded using *MMEMemory:LOAD:STATE* .

Parameters **filename** – is the name of the settings file without path and extension

local_display(*state: bool*) → None

Enable or disable local display of the scope.

Parameters **state** – is the desired local display state

prepare_ultra_segmentation() → None

Make ready for a new acquisition in ultra segmentation mode. This function does one acquisition without ultra segmentation to clear the history and prepare for a new measurement.

read_measurement(*meas_n: int, name: str*) → float

Parameters

- **meas_n** – measurement number 1..8
- **name** – measurement name, for example “MAX”

Returns measured value

run_continuous_acquisition() → None

Start acquiring continuously.

run_single_acquisition() → None

Start a single or Nx acquisition.

save_configuration(*filename: str*) → None

Save the current oscilloscope settings to a file. The filename has to be specified without path and ‘.dfl’ extension, the file will be saved to the configured settings directory.

Information from the manual *SAVe* stores the current instrument settings under the specified number in an intermediate memory. The settings can be recalled using the command **RCL* with the associated number. To transfer the stored instrument settings to a file, use *MMEMemory:STORE:STATE* .

Parameters **filename** – is the name of the settings file without path and extension

save_waveform_history(*filename: str, channel: int, waveform: int = 1*) → None

Save the history of one channel and one waveform to a .bin file. This function is used after an acquisition using sequence trigger mode (with or without ultra segmentation) was performed.

Parameters

- **filename** – is the name (without extension) of the file
- **channel** – is the channel number
- **waveform** – is the waveform number (typically 1)

Raises `RT01024Error` – if storing waveform times out

set_acquire_length(*timerange*: float) → None

Defines the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 … 500 [s]
- Increment: 1E-12 [s]
- *RST = 0.5 [s]

Parameters **timerange** – is the time for one acquisition. Range: 250e-12 … 500 [s]

set_channel_offset(*channel*: int, *offset*: float) → None

Sets the voltage offset of the indicated channel.

- Range: Dependent on the channel scale and coupling [V]
- Increment: Minimum 0.001 [V], may be higher depending on the channel scale and coupling
- *RST = 0

Parameters

- **channel** – is the channel number (1..4)
- **offset** – Offset voltage. Positive values move the waveform down, negative values move it up.

set_channel_position(*channel*: int, *position*: float) → None

Sets the vertical position of the indicated channel as a graphical value.

- Range: -5.0 … 5.0 [div]
- Increment: 0.02
- *RST = 0

Parameters

- **channel** – is the channel number (1..4)
- **position** – is the position. Positive values move the waveform up, negative values move it down.

set_channel_range(*channel*: int, *v_range*: float) → None

Sets the voltage range across the 10 vertical divisions of the diagram. Use the command alternatively instead of `set_channel_scale`.

- Range for range: Depends on attenuation factors and coupling. With 1:1 probe and external attenuations and 50 input coupling, the range is 10 mV to 10 V. For 1 M input coupling, it is 10 mV to 100 V. If the probe and/or external attenuation is changed, multiply the range values by the attenuation factors.
- Increment: 0.01
- *RST = 0.5

Parameters

- **channel** – is the channel number (1..4)
- **v_range** – is the vertical range [V]

set_channel_scale(*channel: int, scale: float*) → None

Sets the vertical scale for the indicated channel. The scale value is given in volts per division.

- Range for scale: depends on attenuation factor and coupling. With 1:1 probe and external attenuations and 50 input coupling, the vertical scale (input sensitivity) is 1 mV/div to 1 V/div. For 1 M input coupling, it is 1 mV/div to 10 V/div. If the probe and/or external attenuation is changed, multiply the values by the attenuation factors to get the actual scale range.
- Increment: 1e-3
- *RST = 0.05

See also: [set_channel_range](#)

Parameters

- **channel** – is the channel number (1..4)
- **scale** – is the vertical scaling [V/div]

set_channel_state(*channel: int, state: bool*) → None

Switches the channel signal on or off.

Parameters

- **channel** – is the input channel (1..4)
- **state** – is True for on, False for off

set_reference_point(*percentage: int*) → None

Sets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]
- *RST = 50 [%]

Parameters **percentage** – is the reference in %

set_repetitions(*number: int*) → None

Set the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- *RST = 1

Parameters **number** – is the number of waveforms to acquire

set_trigger_level(*channel: int, level: float, event_type: int = 1*) → None

Sets the trigger level for the specified event and source.

- Range: -10 to 10 V
- Increment: 1e-3 V
- *RST = 0 V

Parameters

- **channel** – indicates the trigger source.

- 1..4 = channel 1 to 4, available for all event types 1..3
 - 5 = external trigger input on the rear panel for analog signals, available for A-event type
= 1
 - 6..9 = not available
- **level** – is the voltage for the trigger level in [V].
 - **event_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

set_trigger_mode(*mode*: Union[str, hvl_ccb.dev.rs_rto1024.RTO1024.TriggerModes]) → None
Sets the trigger mode which determines the behavior of the instrument if no trigger occurs.

Parameters **mode** – is either auto, normal, or freerun.

Raises **RTO1024Error** – if an invalid triggermode is selected

set_trigger_source(*channel*: int, *event_type*: int = 1) → None
Set the trigger (Event A) source channel.

Parameters

- **channel** – is the channel number (1..4)
- **event_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

start() → None
Start the RTO1024 oscilloscope and bring it into a defined state and remote mode.

stop() → None
Stop the RTO1024 oscilloscope, reset events and close communication. Brings back the device to a state where local operation is possible.

stop_acquisition() → None
Stop any acquisition.

class RT01024Config(*waveforms_path*: str, *settings_path*: str, *backup_path*: str, *spoll_interval*: Union[int, float] = 0.5, *spoll_start_delay*: Union[int, float] = 2, *command_timeout_seconds*: Union[int, float] = 60, *wait_sec_short_pause*: Union[int, float] = 0.1, *wait_sec_enable_history*: Union[int, float] = 1, *wait_sec_post_acquisition_start*: Union[int, float] = 2)
Bases: **hvl_ccb.dev.visa.VisaDeviceConfig**, **hvl_ccb.dev.rs_rto1024._RT01024ConfigDefaultsBase**, **hvl_ccb.dev.rs_rto1024._RT01024ConfigBase**

Configdataclass for the RTO1024 device.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
 Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception RT01024Error

Bases: Exception

class RT01024VisaCommunication(configuration)

Bases: [hvl_ccb.comm.visa.VisaCommunication](#)

Specialization of VisaCommunication for the RTO1024 oscilloscope

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

**class RT01024VisaCommunicationConfig(host: str, interface_type: Union[str,
 hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType] =
 InterfaceType.TCPPIP_INSTR, board: int = 0, port: int = 5025,
 timeout: int = 5000, chunk_size: int = 204800, open_timeout: int =
 1000, write_termination: str = '\n', read_termination: str = '\n',
 visa_backend: str = '')**

Bases: [hvl_ccb.comm.visa.VisaCommunicationConfig](#)

Configuration dataclass for VisaCommunication with specifications for the RTO1024 device class.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

interface_type: Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType] = 2

Interface type of the VISA connection, being one of InterfaceType.

classmethod keys()

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults()

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.


Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

hvl_ccb.dev.se_il2t

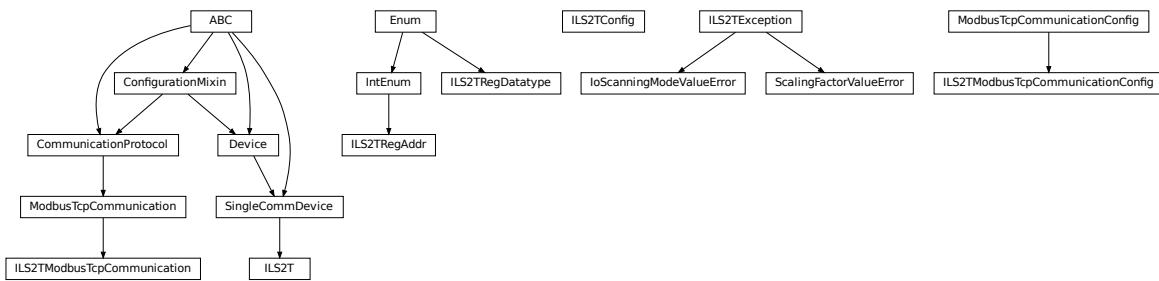

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

class ILS2T(com, dev_config=None)

Bases: [hvl_ccb.dev.base.SingleCommDevice](#)

Schneider Electric ILS2T stepper drive class.

ACTION_JOG_VALUE = 0

The single action value for *ILS2T.Mode.JOG*

class ActionsPtp(value)

Bases: [enum.IntEnum](#)

Allowed actions in the point to point mode (*ILS2T.Mode.PTP*).

ABSOLUTE_POSITION = 0

RELATIVE_POSITION_MOTOR = 2

RELATIVE_POSITION_TARGET = 1

DEFAULT_IO_SCANNING_CONTROL_VALUES = { 'action': 2, 'continue_after_stop_cu': 0, 'disable_driver_di': 0, 'enable_driver_en': 0, 'execute_stop_sh': 0, 'fault_reset_fr': 0, 'mode': 3, 'quick_stop_qs': 0, 'ref_16': 1500, 'ref_32': 0, 'reset_stop_ch': 0 }

Default IO Scanning control mode values

class Mode(value)

Bases: [enum.IntEnum](#)

ILS2T device modes

JOG = 1

PTP = 3

class Ref16Jog(value)

Bases: [enum.Flag](#)

Allowed values for ILS2T ref_16 register (the shown values are the integer representation of the bits), all in Jog mode = 1

```

FAST = 4
NEG = 2
NEG_FAST = 6
NONE = 0
POS = 1
POS_FAST = 5

RegAddr
    Modbus Register Adresses
    alias of hvl\_ccb.dev.se\_ilst2t.ILS2TRegAddr

RegDatatype
    Modbus Register Datatypes
    alias of hvl\_ccb.dev.se\_ilst2t.ILS2TRegDatatype

class State(value)
    Bases: enum.IntEnum
    State machine status values
    ON = 6
    QUICKSTOP = 7
    READY = 4

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

disable(log_warn: bool = True, wait_sec_max: Optional[int] = None) → bool
    Disable the driver of the stepper motor and enable the brake.

    Note: the driver cannot be disabled if the motor is still running.

Parameters

- log_warn – if log a warning in case the motor cannot be disabled.
- wait_sec_max – maximal wait time for the motor to stop running and to disable it; by default, with None, use a config value

Returns True if disable request could and was sent, False otherwise.

do_ioscanning_write(**kwargs: int) → None
    Perform a write operation using IO Scanning mode.

    Parameters kwargs – Keyword-argument list with options to send, remaining are taken from the defaults.

enable() → None
    Enable the driver of the stepper motor and disable the brake.

```

execute_absolute_position(*position: int*) → bool

Execute a absolute position change, i.e. enable motor, perform absolute position change, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

Parameters **position** – absolute position of motor in user defined steps.

Returns *True* if actual position is as expected, *False* otherwise.

execute_relative_step(*steps: int*) → bool

Execute a relative step, i.e. enable motor, perform relative steps, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

Parameters **steps** – Number of steps.

Returns *True* if actual position is as expected, *False* otherwise.

get_dc_volt() → float

Read the DC supply voltage of the motor.

Returns DC input voltage.

get_error_code() → Dict[int, Dict[str, Any]]

Read all messages in fault memory. Will read the full error message and return the decoded values. At the end the fault memory of the motor will be deleted. In addition, *reset_error* is called to re-enable the motor for operation.

Returns Dictionary with all information

get_position() → int

Read the position of the drive and store into status.

Returns Position step value

get_status() → Dict[str, int]

Perform an IO Scanning read and return the status of the motor.

Returns dict with status information.

get_temperature() → int

Read the temperature of the motor.

Returns Temperature in degrees Celsius.

jog_run(*direction: bool = True, fast: bool = False*) → None

Slowly turn the motor in positive direction.

jog_stop() → None

Stop turning the motor in Jog mode.

quickstop() → None

Stops the motor with high deceleration rate and falls into error state. Reset with *reset_error* to recover into normal state.

reset_error() → None

Resets the motor into normal state after quick stop or another error occurred.

set_jog_speed(*slow: int = 60, fast: int = 180*) → None

Set the speed for jog mode. Default values correspond to startup values of the motor.

Parameters

- **slow** – RPM for slow jog mode.

- **fast** – RPM for fast jog mode.

set_max_acceleration(*rpm_minute*: int) → None

Set the maximum acceleration of the motor.

Parameters **rpm_minute** – revolution per minute per minute

set_max_deceleration(*rpm_minute*: int) → None

Set the maximum deceleration of the motor.

Parameters **rpm_minute** – revolution per minute per minute

set_max_rpm(*rpm*: int) → None

Set the maximum RPM.

Parameters **rpm** – revolution per minute (0 < rpm <= RPM_MAX)

Raises **ILS2TException** – if RPM is out of range

set_ramp_type(*ramp_type*: int = -1) → None

Set the ramp type. There are two options available: 0: linear ramp -1: motor optimized ramp

Parameters **ramp_type** – 0: linear ramp | -1: motor optimized ramp

start() → None

Start this device.

stop() → None

Stop this device. Disables the motor (applies brake), disables access and closes the communication protocol.

user_steps(*steps*: int = 16384, *revolutions*: int = 1) → None

Define steps per revolution. Default is 16384 steps per revolution. Maximum precision is 32768 steps per revolution.

Parameters

- **steps** – number of steps in *revolutions*.
- **revolutions** – number of revolutions corresponding to *steps*.

write_absolute_position(*position*: int) → None

Write instruction to turn the motor until it reaches the absolute position. This function does not enable or disable the motor automatically.

Parameters **position** – absolute position of motor in user defined steps.

write_relative_step(*steps*: int) → None

Write instruction to turn the motor the relative amount of steps. This function does not enable or disable the motor automatically.

Parameters **steps** – Number of steps to turn the motor.

```
class ILS2TConfig(rpm_max_init: numbers.Integral = 1500, wait_sec_post_enable: Union[int, float] = 1,
                   wait_sec_max_disable: Union[int, float] = 10, wait_sec_post_cannot_disable: Union[int, float] = 1,
                   wait_sec_post_relative_step: Union[int, float] = 2,
                   wait_sec_post_absolute_position: Union[int, float] = 2)
```

Bases: object

Configuration for the ILS2T stepper motor device.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

rpm_max_init: numbers.Integral = 1500

initial maximum RPM for the motor, can be set up to 3000 RPM. The user is allowed to set a new max RPM at runtime using *ILS2T.set_max_rpm()*, but the value must never exceed this configuration setting.

wait_sec_max_disable: Union[int, float] = 10**wait_sec_post_absolute_position: Union[int, float] = 2****wait_sec_post_CANNOT_DISABLE: Union[int, float] = 1****wait_sec_post_enable: Union[int, float] = 1****wait_sec_post_relative_step: Union[int, float] = 2****exception ILS2TException**

Bases: *Exception*

Exception to indicate problems with the SE ILS2T stepper motor.

class ILS2TModbusTcpCommunication(*configuration*)

Bases: *hvl_ccb.comm.modbus_tcp.ModbusTcpCommunication*

Specific implementation of Modbus/TCP for the Schneider Electric ILS2T stepper motor.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class ILS2TModbusTcpCommunicationConfig(*host: str, unit: int = 255, port: int = 502*)

Bases: *hvl_ccb.comm.modbus_tcp.ModbusTcpCommunicationConfig*

Configuration dataclass for Modbus/TCP communication specific for the Schneider Electric ILS2T stepper motor.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

unit: int = 255

The unit has to be 255 such that IO scanning mode works.

class ILS2TRegAddr(*value*)

Bases: enum.IntEnum

Modbus Register Adresses for Schneider Electric ILS2T stepper drive.

ACCESS_ENABLE = 282

FLT_INFO = 15362

FLT_MEM_DEL = 15112

FLT_MEM_RESET = 15114

IO_SCANNING = 6922

JOGN_FAST = 10506

JOGN_SLOW = 10504

POSITION = 7706

RAMP_ACC = 1556

RAMP_DECEL = 1558

RAMP_N_MAX = 1554

RAMP_TYPE = 1574

SCALE = 1550

TEMP = 7200

VOLT = 7198

```
class ILS2TRegDatatype(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
```

Bases: `aenum.Enum`

Modbus Register Datatypes for Schneider Electric ILS2T stepper drive.

From the manual of the drive:

datatype	byte	min	max
INT8	1 Byte	-128	127
UINT8	1 Byte	0	255
INT16	2 Byte	-32_768	32_767
UINT16	2 Byte	0	65_535
INT32	4 Byte	-2_147_483_648	2_147_483_647
UINT32	4 Byte	0	4_294_967_295
BITS	just 32bits	N/A	N/A

```
INT32 = (-2147483648, 2147483647)
```

```
is_in_range(value: int) → bool
```

```
exception IoScanningModeValueError
```

Bases: `hvl_ccb.dev.se_ilst2t.ILS2TException`


Exception to indicate that the selected IO scanning mode is invalid.

```
exception ScalingFactorValueError
```

Bases: `hvl_ccb.dev.se_ilst2t.ILS2TException`

Exception to indicate that a scaling factor value is invalid.

hvl_ccb.dev.sst_luminox

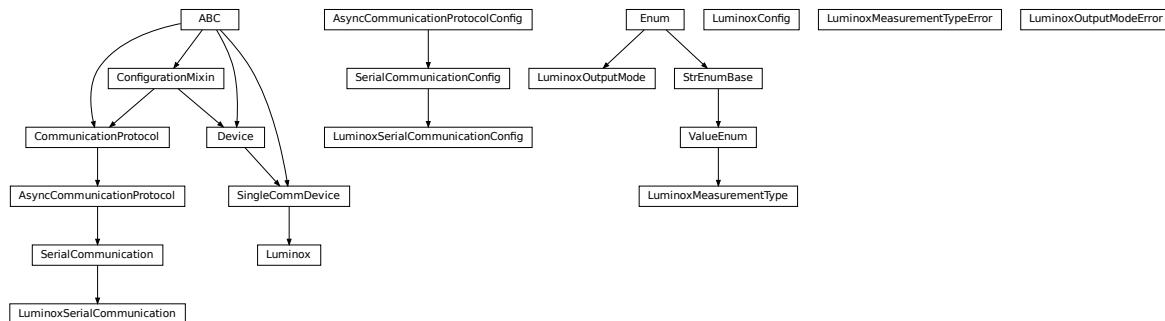

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.

Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

class Luminox(*com, dev_config=None*)
Bases: *hvl_ccb.dev.base.SingleCommDevice*

Luminox oxygen sensor device class.

activate_output(*mode: hvl_ccb.dev.sst_luminox.LuminoxOutputMode*) → None
activate the selected output mode of the Luminox Sensor. :param mode: polling or streaming

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

query_polling(*measurement: Union[str, hvl_ccb.dev.sst_luminox.LuminoxMeasurementType]*) → Union[Dict[Union[str, *hvl_ccb.dev.sst_luminox.LuminoxMeasurementType*], Union[float, int, str]], float, int, str]
Query a value or values of Luminox measurements in the polling mode, according to a given measurement type.

Parameters **measurement** – type of measurement

Returns value of requested measurement

Raises

- **ValueError** – when a wrong key for LuminoxMeasurementType is provided
- **LuminoxOutputModeError** – when polling mode is not activated
- **LuminoxMeasurementTypeError** – when expected measurement value is not read

read_streaming() → Dict[Union[str, *hvl_ccb.dev.sst_luminox.LuminoxMeasurementType*], Union[float, int, str]]
Read values of Luminox in the streaming mode. Convert the single string into separate values.

Returns dictionary with *LuminoxMeasurementType.all_measurements_types()* keys and accordingly type-parsed values.

Raises

- **LuminoxOutputModeError** – when streaming mode is not activated
- **LuminoxMeasurementTypeError** – when any of expected measurement values is not read

start() → None
Start this device. Opens the communication protocol.

stop() → None
Stop the device. Closes also the communication protocol.

class LuminoxConfig(*wait_sec_post_activate: Union[int, float] = 0.5, wait_sec_trials_activate: Union[int, float] = 0.1, nr_trials_activate: int = 5*)
Bases: *object*

Configuration for the SST Luminox oxygen sensor.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys()** → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

nr_trials_activate: int = 5**classmethod optional_defaults()** → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

wait_sec_post_activate: Union[int, float] = 0.5**wait_sec_trials_activate: Union[int, float] = 0.1**

class LuminoxMeasurementType(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*,
boundary=None)

Bases: *hvl_ccb.utils.enum.ValueEnum*

Measurement types for *LuminoxOutputMode.polling*.

The *all_measurements* type will read values for the actual measurement types as given in *LuminoxOutputMode.all_measurements_types()*; it parses multiple single values using regexp's for other measurement types, therefore, no regexp is defined for this measurement type.

all_measurements = 'A'**classmethod all_measurements_types()** → Tuple[*hvl_ccb.dev.sst_luminox.LuminoxMeasurementType*, ...]

A tuple of *LuminoxMeasurementType* enum instances which are actual measurements, i.e. not date of manufacture or software revision.

barometric_pressure = 'P'**property command: str****date_of_manufacture = '# 0'****parse_read_measurement_value**(*read_txt: str*) → Union[Dict[Union[str, *hvl_ccb.dev.sst_luminox.LuminoxMeasurementType*], Union[float, int, str]], float, int, str]**partial_pressure_o2 = '0'**

```
percent_o2 = '%'
sensor_status = 'e'
serial_number = '# 1'
software_revision = '# 2'
temperature_sensor = 'T'
```

LuminoxMeasurementTypeDict

A typing hint for a dictionary holding LuminoxMeasurementType values. Keys are allowed as strings because *LuminoxMeasurementType* is of a *StrEnumBase* type.

alias of Dict[Union[str, *LuminoxMeasurementType*], Union[float, int, str]]

exception LuminoxMeasurementTypeError

Bases: Exception

Wrong measurement type for requested data

LuminoxMeasurementTypeValue

A typing hint for all possible LuminoxMeasurementType values as read in either streaming mode or in a polling mode with *LuminoxMeasurementType.all_measurements*.

Beware: has to be manually kept in sync with *LuminoxMeasurementType* instances *cast_type* attribute values.

alias of Union[float, int, str]

class LuminoxOutputMode(*value*)

Bases: enum.Enum

output mode.

polling = 1

streaming = 0

exception LuminoxOutputModeError

Bases: Exception

Wrong output mode for requested data

class LuminoxSerialCommunication(*configuration*)

Bases: *hvl_ccb.comm.serial.SerialCommunication*

Specific communication protocol implementation for the SST Luminox oxygen sensor. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class LuminoxSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, port:
                                         Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                         Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
                                         <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                         <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                         <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                         Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for SST Luminox is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

SST Luminox does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1

SST Luminox does use one stop bit


terminator: bytes = b'\r\n'

The terminator is CR LF

timeout: Union[int, float] = 3

use 3 seconds timeout as default

hvl_ccb.dev.technix

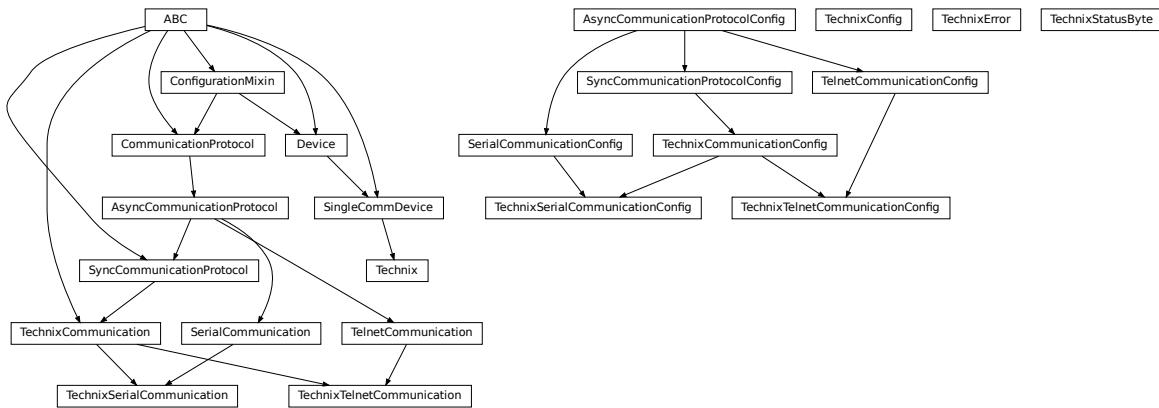

Device classes for “RS 232” and “Ethernet” Interfaces which are used to control power supplies from Technix. Manufacturer homepage: <https://www.technix-hv.com>

The Regulated power Supplies Series and Capacitor Chargers Series from Technix are series of low and high voltage direct current power supplies as well as capacitor chargers. The class *Technix* is tested with a CCR10KV-7,5KJ via an ethernet connection as well as a CCR15-P-2500-OP via a serial connection. Check the code carefully before using it with other devices or device series

This Python package may support the following interfaces from Technix:

- Remote Interface RS232
- Ethernet Remote Interface
- Optic Fiber Remote Interface

```

class Technix(com, dev_config)
    Bases: hvl_ccb.dev.base.SingleCommDevice

    static config_cls()
        Return the default configdataclass class.

        Returns a reference to the default configdataclass class

    property current: Union[int, float]

    default_com_cls() → Union[Type[hvl_ccb.dev.technix.TechinxSerialCommunication],
                                Type[hvl_ccb.dev.technix.TechinxTelnetCommunication]]
        Get the class for the default communication protocol used with this device.

        Returns the type of the standard communication protocol for this device

    get_status_byte() → hvl_ccb.dev.technix.TechinxStatusByte

    property hv: Optional[bool]
    property inhibit: Optional[bool]
    property max_current: Union[int, float]
    property max_voltage: Union[int, float]
    property remote: Optional[bool]
  
```

```
start()
    Open the associated communication protocol.

stop()
    Close the associated communication protocol.

property voltage: Union[int, float]
property voltage_regulation: Optional[bool]

class TechnixCommunication(config)
    Bases: hvl_ccb.comm.base.SyncCommunicationProtocol1, abc.ABC
    Generic communication class for Technix, which can be implemented via TechnixSerialCommunication or TechnixTelnetCommunication

    query(command: str) → str
        Send a command to the interface and handle the status message. Eventually raises an exception.

            Parameters command – Command to send
            Raises TechnixError – if the connection is broken
            Returns Answer from the interface

    class TechnixCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                    encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                    Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int =
                                    10)
        Bases: hvl_ccb.comm.base.SyncCommunicationProtocolConfig

        force_value(fieldname, value)
            Forces a value to a dataclass field despite the class being frozen.

            NOTE: you can define post_force_value method with same signature as this method to do extra processing
            after value has been forced on fieldname.

            Parameters
                • fieldname – name of the field
                • value – value to assign

        classmethod keys() → Sequence[str]
            Returns a list of all configdataclass fields key-names.

            Returns a list of strings containing all keys.

        classmethod optional_defaults() → Dict[str, object]
            Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
            on instantiation.

            Returns a list of strings containing all optional keys.

        classmethod required_keys() → Sequence[str]
            Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
            instantiation.

            Returns a list of strings containing all required keys.

terminator: bytes = b'\r'
    The terminator is CR
```

```
class TechnixConfig(communication_channel: Union[Type[hvl_ccb.dev.technix.TechnixSerialCommunication], Type[hvl_ccb.dev.technix.TechnixTelnetCommunication]], max_voltage: Union[int, float], max_current: Union[int, float], polling_interval_sec: Union[int, float] = 4, post_stop_pause_sec: Union[int, float] = 1, register_pulse_time: Union[int, float] = 0.1)
```

Bases: object

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

```
communication_channel: Union[Type[hvl_ccb.dev.technix.TechnixSerialCommunication], Type[hvl_ccb.dev.technix.TechnixTelnetCommunication]]
```

communication channel between computer and Technix

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

max_current: Union[int, float]

Maximal Output current

max_voltage: Union[int, float]

Maximal Output voltage

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

polling_interval_sec: Union[int, float] = 4

Polling interval in s to maintain to watchdog of the device

post_stop_pause_sec: Union[int, float] = 1

Time to wait after stopping the device

register_pulse_time: Union[int, float] = 0.1

Time for pulsing a register

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception TechnixError

Bases: Exception

Technix related errors.

```
class TechnixSerialCommunication(configuration)
    Bases:           hvl_ccb.dev.technix.TechnixCommunication,      hvl_ccb.comm.serial.
                  SerialCommunication

    static config_cls()
        Return the default configdataclass class.

        Returns a reference to the default configdataclass class

class TechnixSerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, port:
                                         Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                         Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
                                         <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
                                         float, hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                         <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                         <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                         Union[int, float] = 2)
    Bases:           hvl_ccb.dev.technix.TechnixCommunicationConfig,      hvl_ccb.comm.serial.
                  SerialCommunicationConfig

    force_value(fieldname, value)
        Forces a value to a dataclass field despite the class being frozen.

        NOTE: you can define post_force_value method with same signature as this method to do extra processing
              after value has been forced on fieldname.

        Parameters
            • fieldname – name of the field
            • value – value to assign

    classmethod keys() → Sequence[str]
        Returns a list of all configdataclass fields key-names.

        Returns a list of strings containing all keys.

    classmethod optional_defaults() → Dict[str, object]
        Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
        on instantiation.

        Returns a list of strings containing all optional keys.

    classmethod required_keys() → Sequence[str]
        Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
        instantiation.

        Returns a list of strings containing all required keys.

class TechnixStatusByte(value: int)
    Bases: object

    msb_first(idx: int) → Optional[bool]
        Give the Bit at position idx with MSB first

        Parameters idx – Position of Bit as 1...8

        Returns
```

```
class TechnixTelnetCommunication(configuration)
Bases: hvl_ccb.comm.telnet.TelnetCommunication, hvl_ccb.dev.technix.TechinxCommunication
```

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class TechnixTelnetCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, host: Union[str, NoneType] = None, port: int = 4660, timeout: Union[int, float] = 0.2)
```

Bases: *hvl_ccb.comm.telnet.TelnetCommunicationConfig*, *hvl_ccb.dev.technix.TechinxCommunicationConfig*

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 4660

Port at which Technix is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

hvl_ccb.dev.utils

Poller

```
class Poller(spoll_handler: Callable, polling_delay_sec: Union[int, float] = 0, polling_interval_sec: Union[int, float] = 1, polling_timeout_sec: Optional[Union[int, float]] = None)
Bases: object
```

Poller class wrapping `concurrent.futures.ThreadPoolExecutor` which enables passing of results and errors out of the polling thread.

is_polling() → bool

Check if device status is being polled.

Returns *True* when polling thread is set and alive

start_polling() → bool

Start polling.

Returns *True* if was not polling before, *False* otherwise

stop_polling() → bool

Stop polling.

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

Returns *True* if was polling before, *False* otherwise, and last result of the polling function call.


Raises polling function exceptions

wait_for_polling_result()

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

Returns polling function result

Raises polling function errors

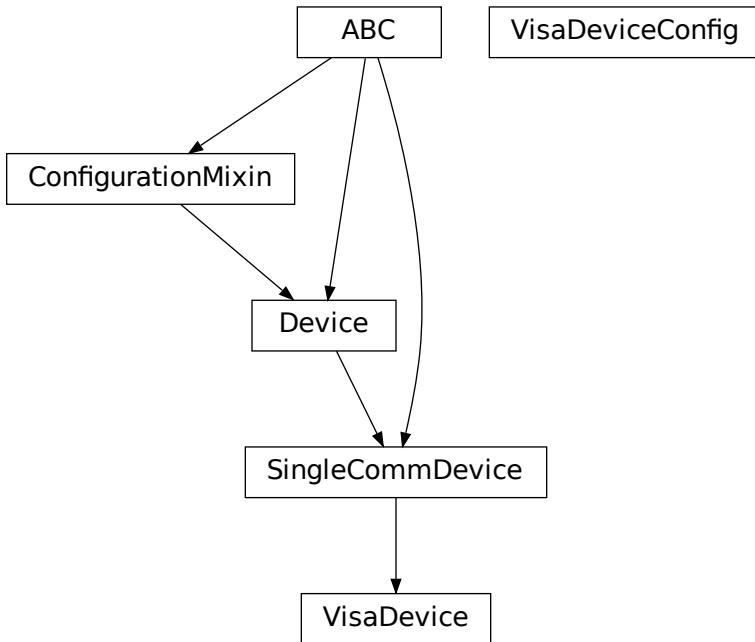hvl_ccb.dev.visa

```
class VisaDevice(com: Union[hvl_ccb.comm.visa.VisaCommunication,
    hvl_ccb.comm.visa.VisaCommunicationConfig, dict], dev_config:
    Optional[Union[hvl_ccb.dev.visa.VisaDeviceConfig, dict]] = None)
Bases: hvl_ccb.dev.base.SingleCommDevice
```

Device communicating over the VISA protocol using VisaCommunication.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls() → Type[hvl_ccb.comm.visa.VisaCommunication]

Return the default communication protocol for this device type, which is VisaCommunication.

Returns the VisaCommunication class

get_error_queue() → str

Read out error queue and logs the error.

Returns Error string

get_identification() → str

Queries “*IDN?” and returns the identification string of the connected device.

Returns the identification string of the connected device

reset() → None

Send “*RST” and “*CLS” to the device. Typically sets a defined state.

spoll_handler()

Reads the status byte and decodes it. The status byte STB is defined in IEEE 488.2. It provides a rough overview of the instrument status.

Returns**start() → None**

Start the VisaDevice. Sets up the status poller and starts it.

Returns**stop() → None**

Stop the VisaDevice. Stops the polling thread and closes the communication protocol.

Returns**wait_operation_complete(timeout: Optional[float] = None) → bool**

Waits for a operation complete event. Returns after timeout [s] has expired or the operation complete event has been caught.

Parameters **timeout** – Time in seconds to wait for the event; *None* for no timeout.

Returns True, if OPC event is caught, False if timeout expired

class VisaDeviceConfig(spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2)
Bases: hvl_ccb.dev.visa._VisaDeviceConfigDefaultsBase, hvl_ccb.dev.visa._VisaDeviceConfigBase

Configdataclass for a VISA device.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.


Module contents

Devices subpackage.

`hvl_ccb.utils`

Submodules

`hvl_ccb.utils.enum`

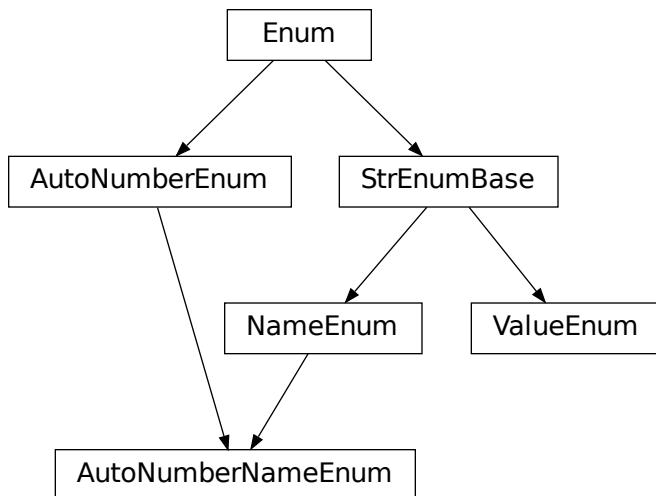

class `AutoNumberNameEnum`(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.NameEnum`, `aenum.AutoNumberEnum`

Auto-numbered enum with names used as string representation, and with lookup and equality based on this representation.

class `NameEnum`(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.StrEnumBase`

Enum with names used as string representation, and with lookup and equality based on this representation.

class `StrEnumBase`(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `aenum.Enum`

String representation-based equality and lookup.

class `ValueEnum`(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: `hvl_ccb.utils.enum.StrEnumBase`

Enum with string representation of values used as string representation, and with lookup and equality based on this representation.

Attention: to avoid errors, best use together with *unique* enum decorator.

hvl_ccb.utils.typing

Additional Python typing module utilities

ConvertibleTypes

Typing hint for data type that can be used in conversion

alias of Union[int, float, List[Union[int, float]], Tuple[Union[int, float]], Dict[str, Union[int, float]], numpy.ndarray]

Number

Typing hint auxiliary for a Python base number types: *int* or *float*.

alias of Union[int, float]

check_generic_type(value, type_, name='instance')

Check if *value* is of a generic type *type_*. Raises *TypeError* if it's not.

Parameters

- **name** – name to report in case of an error
- **value** – value to check
- **type** – generic type to check against

is_generic_type_hint(type_)

Check if class is a generic type, for example *Union[int, float]*, *List[int]*, or *List*.

Parameters **type** – type to check

hvl_ccb.utils.validation

validate_bool(x_name: str, x: object, logger: Optional[logging.Logger] = None) → None

Validate if given input *x* is a *bool*.

Parameters

- **x_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as boolean
- **logger** – logger of the calling submodule

Raises **TypeError** – when the validated input does not have boolean type

validate_number(x_name: str, x: object, limits: Optional[Tuple] = (None, None), number_type:

Union[Type[Union[int, float]], Tuple[Type[Union[int, float]], ...]] = (<class 'int'>, <class 'float'>), logger: Optional[logging.Logger] = None) → None

Validate if given input *x* is a number of given *number_type* type, with value between given *limits[0]* and *limits[1]* (inclusive), if not *None*.

Parameters

- **x_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as number of given type within given range
- **logger** – logger of the calling submodule
- **limits** – [lower, upper] limit, with *None* denoting no limit: [-inf, +inf]

- **number_type** – expected type or tuple of types of a number, by default (*int, float*)


Raises

- **TypeError** – when the validated input does not have expected type
- **ValueError** – when the validated input has correct number type but is not within given range

Module contents

4.1.2 Submodules

hvl_ccb.configuration

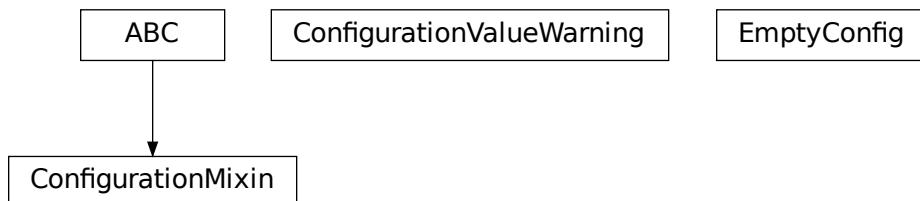

providing classes for handling configuration for communication protocols and devices.

Facilities

class ConfigurationMixin(configuration)

Bases: abc.ABC

Mixin providing configuration to a class.

property config

ConfigDataclass property.

Returns the configuration

abstract static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

configuration_save_json(path: str) → None

Save current configuration as JSON file.

Parameters **path** – path to the JSON file.

classmethod from_json(filename: str)

Instantiate communication protocol using configuration from a JSON file.

Parameters **filename** – Path and filename to the JSON configuration

exception ConfigurationValueWarning

Bases: UserWarning

User warnings category for values of @configdataclass fields.

class EmptyConfig

Bases: object

Empty configuration dataclass.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns a list of strings containing all required keys.

configdataclass(direct_decoration=None, frozen=True)

Decorator to make a class a configdataclass. Types in these dataclasses are enforced. Implement a function *clean_values(self)* to do additional checking on value ranges etc.

It is possible to inherit from a configdataclass and re-decorate it with @configdataclass. In a subclass, default values can be added to existing fields. Note: adding additional non-default fields is prone to errors, since the order has to be respected through the whole chain (first non-default fields, only then default-fields).

Parameters **frozen** – defaults to True. False allows to later change configuration values. Attention:
if configdataclass is not frozen and a value is changed, typing is not enforced anymore!
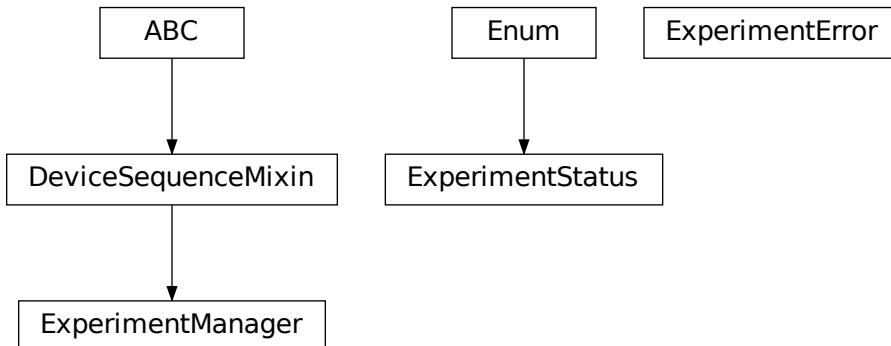
`hvl_ccb.experiment_manager`

Main module containing the top level `ExperimentManager` class. Inherit from this class to implement your own experiment functionality in another project and it will help you start, stop and manage your devices.

`exception ExperimentError`

Bases: `Exception`

Exception to indicate that the current status of the experiment manager is on `ERROR` and thus no operations can be made until reset.

`class ExperimentManager(*args, **kwargs)`
Bases: `hvl_ccb.dev.base.DeviceSequenceMixin`

Experiment Manager can start and stop communication protocols and devices. It provides methods to queue commands to devices and collect results.

`add_device(name: str, device: hvl_ccb.dev.base.Device) → None`

Add a new device to the manager. If the experiment is running, automatically start the device. If a device with this name already exists, raise an exception.

Parameters

- **`name`** – is the name of the device.
- **`device`** – is the instantiated `Device` object.

`Raises DeviceExistingException –`

`devices_failed_start: Dict[str, hvl_ccb.dev.base.Device]`

Dictionary of named device instances from the sequence for which the most recent `start()` attempt failed.

Empty if `stop()` was called last; cf. `devices_failed_stop`.

`devices_failed_stop: Dict[str, hvl_ccb.dev.base.Device]`

Dictionary of named device instances from the sequence for which the most recent `stop()` attempt failed.

Empty if `start()` was called last; cf. `devices_failed_start`.

`finish() → None`

Stop experimental setup, stop all devices.

```
is_error() → bool
    Returns true, if the status of the experiment manager is error.
    Returns True if on error, false otherwise

is_finished() → bool
    Returns true, if the status of the experiment manager is finished.
    Returns True if finished, false otherwise

is_running() → bool
    Returns true, if the status of the experiment manager is running.
    Returns True if running, false otherwise

run() → None
    Start experimental setup, start all devices.

start() → None
    Alias for ExperimentManager.run()

property status: hvl_ccb.experiment_manager.ExperimentStatus
    Get experiment status.
    Returns experiment status enum code.

stop() → None
    Alias for ExperimentManager.finish()

class ExperimentStatus(value)
    Bases: enum.Enum
        Enumeration for the experiment status
        ERROR = 5
        FINISHED = 4
        FINISHING = 3
        INITIALIZED = 0
        INITIALIZING = -1
        RUNNING = 2
        STARTING = 1
```

4.1.3 Module contents

Top-level package for HVL Common Code Base.

CONTRIBUTING

Contributions are welcome, and they are greatly appreciated! Every little bit helps, and credit will always be given.

You can contribute in many ways:

5.1 Types of Contributions

5.1.1 Report Bugs

Report bugs at https://gitlab.com/ethz_hvl/hvl_ccb/issues.

If you are reporting a bug, please include:

- Your operating system name and version.
- Any details about your local setup that might be helpful in troubleshooting.
- Detailed steps to reproduce the bug.

5.1.2 Fix Bugs

Look through the GitLab issues for bugs. Anything tagged with “bug” and “help wanted” is open to whoever wants to implement it.

5.1.3 Implement Features

Look through the GitLab issues for features. Anything tagged with “enhancement” and “help wanted” is open to whoever wants to implement it.

5.1.4 Write Documentation

HVL Common Code Base could always use more documentation, whether as part of the official HVL Common Code Base docs, in docstrings, or even on the web in blog posts, articles, and such.

5.1.5 Submit Feedback

The best way to send feedback is to file an issue at https://gitlab.com/ethz_hvl/hvl_ccb/issues.

If you are proposing a feature:

- Explain in detail how it would work.
- Keep the scope as narrow as possible, to make it easier to implement.
- Remember that this is a volunteer-driven project, and that contributions are welcome :)

5.2 Get Started!

Ready to contribute? Here's how to set up *hvl_ccb* for local development.

1. Clone *hvl_ccb* repo from GitLab.

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

2. Install your local copy into a virtualenv. Assuming you have `virtualenvwrapper` installed, this is how you set up your fork for local development:

```
$ mkvirtualenv hvl_ccb
$ cd hvl_ccb/
$ pip install -e .[all]
$ pip install -r requirements_dev.txt
```

3. Create a branch for local development:

```
$ git checkout -b name-of-your-bugfix-or-feature
```

Now you can make your changes locally.

4. When you're done making changes, check that your changes pass flake8 and the tests, including testing other Python versions with tox:

```
$ flake8 hvl_ccb tests
$ python setup.py test or py.test
$ tox
```

To get flake8 and tox, just pip install them into your virtualenv. You can also use the provided make-like shell script to run flake8 and tests:

```
$ ./make.sh lint
$ ./make.sh test
```

5. Commit your changes and push your branch to GitLab:

```
$ git add .
$ git commit -m "Your detailed description of your changes."
$ git push origin name-of-your-bugfix-or-feature
```

6. Submit a merge request through the GitLab website.

5.3 Merge Request Guidelines

Before you submit a merge request, check that it meets these guidelines:

1. The merge request should include tests.
2. If the merge request adds functionality, the docs should be updated. Put your new functionality into a function with a docstring, and add the feature to the list in README.rst.
3. The merge request should work for Python 3.7. Check https://gitlab.com/ethz_hvl/hvl_ccb/merge_requests and make sure that the tests pass for all supported Python versions.

5.4 Tips

- To run tests from a single file:

```
$ py.test tests/test_hvl_ccb.py
```

or a single test function:

```
$ py.test tests/test_hvl_ccb.py::test_command_line_interface
```

- If your tests are slow, profile them using the pytest-profiling plugin:

```
$ py.test tests/test_hvl_ccb.py --profile
```

or for a graphical overview (you need a SVG image viewer):

```
$ py.test tests/test_hvl_ccb.py --profile-svg
$ open prof/combined.svg
```

- To add dependency, edit appropriate *requirements variable in the setup.py file and re-run:

```
$ python setup.py develop
```

- To generate a PDF version of the Sphinx documentation instead of HTML use:

```
$ rm -rf docs/hvl_ccb.rst docs/modules.rst docs/_build && sphinx-apidoc -o docs/hvl_ccb && python -msphinx -M latexpdf docs/ docs/_build
```

This command can also be run through the make-like shell script:

```
$ ./make.sh docs-pdf
```

This requires a local installation of a LaTeX distribution, e.g. MikTeX.

5.5 Deploying

A reminder for the maintainers on how to deploy. Create `release-N.M.K` branch. Make sure all your changes are committed. Update or create entry in `HISTORY.rst` file, and, if applicable, update `AUTHORS.rst` file, update features tables in `README.rst` file, and update API docs:

```
$ make docs
```

Commit all of the above, except for the `docs/hvl_ccb.dev.picotech_pt104.rst`, and then run:

```
$ bumpversion patch # possible: major / minor / patch
$ git push
$ git push --tags
$ make release
```

Merge the release branch into master and devel branches with `--no-ff` flag and delete the release branch:

```
$ git switch master
$ git merge --no-ff release-N.M.K
$ git push
$ git switch devel
$ git merge --no-ff release-N.M.K
$ git push
$ git push --delete origin release-N.M.K
$ git branch --delete release-N.M.K
```

Finally, go to https://gitlab.com/ethz_hvl/hvl_ccb/tags/, select the latest release tag, press “Edit release notes” and add release notes (corresponding entry from `HISTORY.rst` file, but consider also additional brief header or synopsis if needed).

CREDITS

6.1 Maintainers

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>

6.2 Authors

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- David Gruber <dev@davidgruber.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Alise Chachereau <chachereau@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- David Taylor <dtaylor@ethz.ch>

6.3 Contributors

- Luca Nembrini <lucane@student.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>
- Raphael Faerber <raphael.faerber@eeh.ee.ethz.ch>
- Ruben Stadler <rstadler@student.ethz.ch>
- Joseph Engelbrecht <engelbrecht@eeh.ee.ethz.ch>
- Hanut Vemulapalli <vemulapalli@eeh.ee.ethz.ch>

HISTORY

7.1 0.8.3 (2021-09-27)

- New data conversion functions in `utils.conversion_sensor` and `utils.conversion_unit` modules. Note: to use these functions you must install `hvl_ccb` with extra requirement, either `hvl_ccb[conversion]` or `hvl_ccb[all]`.
- Improved documentation with respect to installation of external libraries.

7.2 0.8.2 (2021-08-27)

- New functionality in `dev.labjack.LabJack`:
 - configure clock and send timed pulse sequences
 - set DAC/analog output voltage
- Bugfix: ignore random bits sent by to `dev.newport.NewportSMC100PP` controller during start-up/powering-up.

7.3 0.8.1 (2021-08-13)

- Add Python version check (min version error; max version warning).
- Daily checks for upstream dependencies compatibility and devel environment improvements.

7.4 0.8.0 (2021-07-02)

- TCP communication protocol.
- Lauda PRO RP 245 E circulation thermostat device over TCP.
- Pico Technology PT-104 Platinum Resistance Data Logger device as a wrapper of the Python bindings for the PicoSDK.
- In `com.visa.VisaCommunication`: periodic status polling when VISA/TCP keep alive connection is not supported by a host.

7.5 0.7.1 (2021-06-04)

- New `utils.validation` submodule with `validate_bool` and `validate_number` utilities extracted from internal use within a `dev.tiepie` subpackage.
- In `comm.serial.SerialCommunication`:
 - strict encoding errors handling strategy for subclasses,
 - user warning for a low communication timeout value.

7.6 0.7.0 (2021-05-25)

- The `dev.tiepie` module was splitted into a subpackage with, in particular, submodules for each of the device types – oscilloscope, generator, and i2c – and with backward-incompatible direct imports from the submodules.
- In `dev.technix`:
 - fixed communication crash on nested status byte query;
 - added enums for GET and SET register commands.
- Further minor logging improvements: added missing module level logger and removed some error logs in `except` blocks used for a flow control.
- In `examples/` folder renamed consistently all the examples.
- In API documentation: fix incorrect links mapping on inheritance diagrams.

7.7 0.6.1 (2021-05-08)

- In `dev.tiepie`:
 - dynamically set oscilloscope's channel limits in `OscilloscopeChannelParameterLimits`: `input_range` and `trigger_level_abs`, incl. update of latter on each change of `input_range` value of a `TiePieOscilloscopeChannelConfig` instances;
 - quick fix for opening of combined instruments by disabling `OscilloscopeParameterLimits`. `trigger_delay` (an advanced feature);
 - enable automatic devices detection to be able to find network devices with `TiePieOscilloscope`. `list_devices()`.
- Fix `examples/example_labjack.py`.
- Improved logging: consistently use module level loggers, and always log exception tracebacks.
- Improve API documentation: separate pages per modules, each with an inheritance diagram as an overview.

7.8 0.6.0 (2021-04-23)

- Technix capacitor charger using either serial connection or Telnet protocol.
- **Extensions, improvements and fixes in existing devices:**
 - In `dev.tiepie.TiePieOscilloscope`:
 - * redesigned measurement start and data collection API, incl. time out argument, with no/infinite time out option;
 - * trigger allows now a no/infinite time out;
 - * record length and trigger level were fixed to accept, respectively, floating point and integer numbers;
 - * fixed resolution validation bug;
 - `dev.heinzinger.HeinzingerDI` and `dev.rs_rto1024.RT01024` instances are now resilient to multiple `stop()` calls.
 - In `dev.crylas.CryLasLaser`: default configuration timeout and polling period were adjusted;
 - Fixed PSI9080 example script.
- **Package and source code improvements:**
 - Update to backward-incompatible `pyvisa-py>=0.5.2`. Developers, do update your local development environments!
 - External libraries, like LibTiePie SDK or LJM Library, are now not installed by default; they are now extra installation options.
 - Added Python 3.9 support.
 - Improved number formatting in logs.
 - Typing improvements and fixes for `mypy>=0.800`.

7.9 0.5.0 (2020-11-11)

- TiePie USB oscilloscope, generator and I2C host devices, as a wrapper of the Python bindings for the LibTiePie SDK.
- a FuG Elektronik Power Supply (e.g. Capacitor Charger HCK) using the built-in ADDAT controller with the Probus V protocol over a serial connection
- All devices polling status or measurements use now a `dev.utils.Poller` utility class.
- **Extensions and improvements in existing devices:**
 - In `dev.rs_rto1024.RT01024`: added Channel state, scale, range, position and offset accessors, and measurements activation and read methods.
 - In `dev.sst_luminox.Luminox`: added querying for all measurements in polling mode, and made output mode activation more robust.
 - In `dev.newport.NewportSMC100PP`: an error-prone `wait_until_move_finished` method of replaced by a fixed waiting time, device operations are now robust to a power supply cut, and device restart is not required to apply a start configuration.
- **Other minor improvements:**

- Single failure-safe starting and stopping of devices sequenced via `dev.base.DeviceSequenceMixin`.
- Moved `read_text_nonempty` up to `comm.serial.SerialCommunication`.
- Added development Dockerfile.
- Updated package and development dependencies: `pymodbus`, `pytest-mock`.

7.10 0.4.0 (2020-07-16)

- **Significantly improved new Supercube device controller:**
 - more robust error-handling,
 - status polling with generic Poller helper,
 - messages and status boards.
 - tested with a physical device,
- Improved OPC UA client wrapper, with better error handling, incl. re-tries on `concurrent.futures.TimeoutError`.
- SST Luminox Oxygen sensor device controller.
- **Backward-incompatible changes:**
 - `CommunicationProtocol.access_lock` has changed type from `threading.Lock` to `threading.RLock`.
 - `ILS2T.relative_step` and `ILS2T.absolute_position` are now called, respectively, `ILS2T.write_relative_step` and `ILS2T.write_absolute_position`.
- **Minor bugfixes and improvements:**
 - fix use of max resolution in `Labjack.set_ain_resolution()`,
 - resolve ILS2T devices relative and absolute position setters race condition,
 - added acoustic horn function in the 2015 Supercube.
- **Toolchain changes:**
 - add Python 3.8 support,
 - drop pytest-runner support,
 - ensure compatibility with `labjack_ljm` 2019 version library.

7.11 0.3.5 (2020-02-18)

- Fix issue with reading integers from LabJack LJM Library (device's product ID, serial number etc.)
- Fix development requirements specification (tox version).

7.12 0.3.4 (2019-12-20)

- New devices using serial connection:
 - Heinzinger Digital Interface I/II and a Heinzinger PNC power supply
 - Q-switched Pulsed Laser and a laser attenuator from CryLas
 - Newport SMC100PP single axis motion controller for 2-phase stepper motors
 - Pfeiffer TPG controller (TPG 25x, TPG 26x and TPG 36x) for Compact pressure Gauges
- PEP 561 compatibility and related corrections for static type checking (now in CI)
- Refactorings:
 - Protected non-thread safe read and write in communication protocols
 - Device sequence mixin: start/stop, add/rm and lookup
 - `.format()` to f-strings
 - more enumerations and a quite some improvements of existing code
- Improved error docstrings (`:raises:` annotations) and extended tests for errors.

7.13 0.3.3 (2019-05-08)

- Use PyPI labjack-ljm (no external dependencies)

7.14 0.3.2 (2019-05-08)

- INSTALLATION.rst with LJMPython prerequisite info

7.15 0.3.1 (2019-05-02)

- readthedocs.org support

7.16 0.3 (2019-05-02)

- Prevent an automatic close of VISA connection when not used.
- Rhode & Schwarz RTO 1024 oscilloscope using VISA interface over `TCP::INSTR`.
- Extended tests incl. messages sent to devices.
- Added Supercube device using an OPC UA client
- Added Supercube 2015 device using an OPC UA client (for interfacing with old system version)

7.17 0.2.1 (2019-04-01)

- Fix issue with LJMPython not being installed automatically with setuptools.

7.18 0.2.0 (2019-03-31)

- LabJack LJM Library communication wrapper and LabJack device.
- Modbus TCP communication protocol.
- Schneider Electric ILS2T stepper motor drive device.
- Elektro-Automatik PSI9000 current source device and VISA communication wrapper.
- Separate configuration classes for communication protocols and devices.
- Simple experiment manager class.

7.19 0.1.0 (2019-02-06)

- Communication protocol base and serial communication implementation.
- Device base and MBW973 implementation.

**CHAPTER
EIGHT**

INDICES AND TABLES

- genindex
- modindex
- search

PYTHON MODULE INDEX

h

hvl_ccb, 186
hvl_ccb.comm, 33
hvl_ccb.comm.base, 11
hvl_ccb.comm.labjack_ljm, 15
hvl_ccb.comm.modbus_tcp, 18
hvl_ccb.comm.opc, 20
hvl_ccb.comm.serial, 23
hvl_ccb.comm.tcp, 26
hvl_ccb.comm.telnet, 28
hvl_ccb.comm.visa, 30
hvl_ccb.configuration, 183
hvl_ccb.dev, 181
hvl_ccb.dev.base, 84
hvl_ccb.dev.crylas, 87
hvl_ccb.dev.ea_psi9000, 96
hvl_ccb.dev.fug, 100
hvl_ccb.dev.heinzinger, 112
hvl_ccb.dev.labjack, 118
hvl_ccb.dev.lauda, 124
hvl_ccb.dev.mbw973, 131
hvl_ccb.dev.newport, 134
hvl_ccb.dev.pfeiffer_tpg, 148
hvl_ccb.dev.rs_rto1024, 154
hvl_ccb.dev.se_ilis2t, 162
hvl_ccb.dev.sst_luminox, 168
hvl_ccb.dev.supercube, 57
hvl_ccb.dev.supercube.base, 33
hvl_ccb.dev.supercube.constants, 39
hvl_ccb.dev.supercube.typ_a, 53
hvl_ccb.dev.supercube.typ_b, 55
hvl_ccb.dev.supercube2015, 72
hvl_ccb.dev.supercube2015.base, 57
hvl_ccb.dev.supercube2015.constants, 62
hvl_ccb.dev.supercube2015.typ_a, 70
hvl_ccb.dev.technix, 173
hvl_ccb.dev.tiepie, 83
hvl_ccb.dev.tiepie.base, 72
hvl_ccb.dev.tiepie.channel, 74
hvl_ccb.dev.tiepie.device, 77
hvl_ccb.dev.tiepie.generator, 78
hvl_ccb.dev.tiepie.i2c, 79

hvl_ccb.dev.tiepie.oscilloscope, 80
hvl_ccb.dev.tiepie.utils, 83
hvl_ccb.dev.utils, 177
hvl_ccb.dev.visa, 179
hvl_ccb.experiment_manager, 185
hvl_ccb.utils, 183
hvl_ccb.utils.enum, 181
hvl_ccb.utils.typing, 182
hvl_ccb.utils.validation, 182

INDEX

A

A (*HeinzingerPNC.UnitCurrent attribute*), 115
A (*SupercubeOpcEndpoint attribute*), 52, 69
ABSOLUTE (*TiePieOscilloscopeTriggerLevelMode attribute*), 76
ABSOLUTE_POSITION (*ILS2T.ActionsPtp attribute*), 162
AC (*NewportConfigCommands attribute*), 134
AC_DoubleStage_150kV (*PowerSetup attribute*), 51, 68
AC_DoubleStage_200kV (*PowerSetup attribute*), 51, 68
AC_SingleStage_100kV (*PowerSetup attribute*), 51, 68
AC_SingleStage_50kV (*PowerSetup attribute*), 51, 68
ACA (*TiePieOscilloscopeChannelCoupling attribute*), 75
acceleration (*NewportSMC100PPConfig attribute*), 142
ACCESS_ENABLE (*ILS2TRegAddr attribute*), 167
access_lock (*CommunicationProtocol attribute*), 14
ACTION_JOG_VALUE (*ILS2T attribute*), 162
activate_measurements() (*RTO1024 method*), 155
activate_output() (*Luminox method*), 169
activated (*BreakdownDetection attribute*), 44, 63
ACTIVE (*CryLasLaser.AnswersStatus attribute*), 90
active (*OpcControl attribute*), 50
actualsetvalue (*FuGProbusVSetRegisters property*), 109
ACV (*TiePieOscilloscopeChannelCoupling attribute*), 75
adc_mode (*FuGProbusVMonitorRegisters property*), 108
add_device() (*DeviceSequenceMixin method*), 85
add_device() (*ExperimentManager method*), 185
ADDR_INCORRECT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 144
address (*NewportSMC100PPConfig attribute*), 142
address (*VisaCommunicationConfig property*), 31
address() (*VisaCommunicationConfig.InterfaceType method*), 31
ADMODE (*FuGProbusIVCommands attribute*), 106
Alarm0 (*AlarmText attribute*), 62
Alarm1 (*Alarms attribute*), 40
Alarm1 (*AlarmText attribute*), 39, 62
Alarm10 (*Alarms attribute*), 40
Alarm10 (*AlarmText attribute*), 39, 62
Alarm100 (*Alarms attribute*), 40

Alarm101 (*Alarms attribute*), 40
Alarm102 (*Alarms attribute*), 40
Alarm103 (*Alarms attribute*), 40
Alarm104 (*Alarms attribute*), 40
Alarm105 (*Alarms attribute*), 40
Alarm106 (*Alarms attribute*), 40
Alarm107 (*Alarms attribute*), 40
Alarm108 (*Alarms attribute*), 40
Alarm109 (*Alarms attribute*), 40
Alarm11 (*Alarms attribute*), 40
Alarm11 (*AlarmText attribute*), 39, 62
Alarm110 (*Alarms attribute*), 41
Alarm111 (*Alarms attribute*), 41
Alarm112 (*Alarms attribute*), 41
Alarm113 (*Alarms attribute*), 41
Alarm114 (*Alarms attribute*), 41
Alarm115 (*Alarms attribute*), 41
Alarm116 (*Alarms attribute*), 41
Alarm117 (*Alarms attribute*), 41
Alarm118 (*Alarms attribute*), 41
Alarm119 (*Alarms attribute*), 41
Alarm12 (*Alarms attribute*), 41
Alarm12 (*AlarmText attribute*), 39, 62
Alarm120 (*Alarms attribute*), 41
Alarm121 (*Alarms attribute*), 41
Alarm122 (*Alarms attribute*), 41
Alarm123 (*Alarms attribute*), 41
Alarm124 (*Alarms attribute*), 41
Alarm125 (*Alarms attribute*), 41
Alarm126 (*Alarms attribute*), 41
Alarm127 (*Alarms attribute*), 41
Alarm128 (*Alarms attribute*), 41
Alarm129 (*Alarms attribute*), 41
Alarm13 (*Alarms attribute*), 41
Alarm13 (*AlarmText attribute*), 39, 62
Alarm130 (*Alarms attribute*), 41
Alarm131 (*Alarms attribute*), 41
Alarm132 (*Alarms attribute*), 41
Alarm133 (*Alarms attribute*), 41
Alarm134 (*Alarms attribute*), 41
Alarm135 (*Alarms attribute*), 41
Alarm136 (*Alarms attribute*), 41

Alarm137 (*Alarms attribute*), 41
Alarm138 (*Alarms attribute*), 41
Alarm139 (*Alarms attribute*), 41
Alarm14 (*Alarms attribute*), 41
Alarm14 (*AlarmText attribute*), 39, 62
Alarm140 (*Alarms attribute*), 41
Alarm141 (*Alarms attribute*), 41
Alarm142 (*Alarms attribute*), 41
Alarm143 (*Alarms attribute*), 42
Alarm144 (*Alarms attribute*), 42
Alarm145 (*Alarms attribute*), 42
Alarm146 (*Alarms attribute*), 42
Alarm147 (*Alarms attribute*), 42
Alarm148 (*Alarms attribute*), 42
Alarm149 (*Alarms attribute*), 42
Alarm15 (*Alarms attribute*), 42
Alarm15 (*AlarmText attribute*), 39
Alarm150 (*Alarms attribute*), 42
Alarm151 (*Alarms attribute*), 42
Alarm16 (*Alarms attribute*), 42
Alarm16 (*AlarmText attribute*), 39
Alarm17 (*Alarms attribute*), 42
Alarm17 (*AlarmText attribute*), 39, 62
Alarm18 (*Alarms attribute*), 42
Alarm18 (*AlarmText attribute*), 39
Alarm19 (*Alarms attribute*), 42
Alarm19 (*AlarmText attribute*), 39, 62
Alarm2 (*Alarms attribute*), 42
Alarm2 (*AlarmText attribute*), 39, 62
Alarm20 (*Alarms attribute*), 42
Alarm20 (*AlarmText attribute*), 39, 62
Alarm21 (*Alarms attribute*), 42
Alarm21 (*AlarmText attribute*), 39, 62
Alarm22 (*Alarms attribute*), 42
Alarm22 (*AlarmText attribute*), 39, 62
Alarm23 (*Alarms attribute*), 42
Alarm23 (*AlarmText attribute*), 39
Alarm24 (*Alarms attribute*), 42
Alarm24 (*AlarmText attribute*), 39
Alarm25 (*Alarms attribute*), 42
Alarm25 (*AlarmText attribute*), 39
Alarm26 (*Alarms attribute*), 42
Alarm26 (*AlarmText attribute*), 40
Alarm27 (*Alarms attribute*), 42
Alarm28 (*Alarms attribute*), 42
Alarm29 (*Alarms attribute*), 42
Alarm3 (*Alarms attribute*), 42
Alarm3 (*AlarmText attribute*), 40, 63
Alarm30 (*Alarms attribute*), 42
Alarm31 (*Alarms attribute*), 42
Alarm32 (*Alarms attribute*), 42
Alarm33 (*Alarms attribute*), 42
Alarm34 (*Alarms attribute*), 42
Alarm35 (*Alarms attribute*), 42
Alarm36 (*Alarms attribute*), 42
Alarm37 (*Alarms attribute*), 42
Alarm38 (*Alarms attribute*), 42
Alarm39 (*Alarms attribute*), 42
Alarm4 (*Alarms attribute*), 43
Alarm4 (*AlarmText attribute*), 40, 63
Alarm40 (*Alarms attribute*), 43
Alarm41 (*Alarms attribute*), 43
Alarm41 (*AlarmText attribute*), 40
Alarm42 (*Alarms attribute*), 43
Alarm42 (*AlarmText attribute*), 40
Alarm43 (*Alarms attribute*), 43
Alarm43 (*AlarmText attribute*), 40
Alarm44 (*Alarms attribute*), 43
Alarm44 (*AlarmText attribute*), 40
Alarm45 (*Alarms attribute*), 43
Alarm45 (*AlarmText attribute*), 40
Alarm46 (*Alarms attribute*), 43
Alarm46 (*AlarmText attribute*), 40
Alarm47 (*Alarms attribute*), 43
Alarm47 (*AlarmText attribute*), 40
Alarm48 (*Alarms attribute*), 43
Alarm48 (*AlarmText attribute*), 40
Alarm49 (*Alarms attribute*), 43
Alarm5 (*Alarms attribute*), 43
Alarm5 (*AlarmText attribute*), 40, 63
Alarm50 (*Alarms attribute*), 43
Alarm51 (*Alarms attribute*), 43
Alarm52 (*Alarms attribute*), 43
Alarm53 (*Alarms attribute*), 43
Alarm54 (*Alarms attribute*), 43
Alarm55 (*Alarms attribute*), 43
Alarm56 (*Alarms attribute*), 43
Alarm57 (*Alarms attribute*), 43
Alarm58 (*Alarms attribute*), 43
Alarm59 (*Alarms attribute*), 43
Alarm6 (*Alarms attribute*), 43
Alarm6 (*AlarmText attribute*), 40, 63
Alarm60 (*Alarms attribute*), 43
Alarm61 (*Alarms attribute*), 43
Alarm62 (*Alarms attribute*), 43
Alarm63 (*Alarms attribute*), 43
Alarm64 (*Alarms attribute*), 43
Alarm65 (*Alarms attribute*), 43
Alarm66 (*Alarms attribute*), 43
Alarm67 (*Alarms attribute*), 43
Alarm68 (*Alarms attribute*), 43
Alarm69 (*Alarms attribute*), 43
Alarm7 (*Alarms attribute*), 43
Alarm7 (*AlarmText attribute*), 40, 63
Alarm70 (*Alarms attribute*), 43
Alarm71 (*Alarms attribute*), 43
Alarm72 (*Alarms attribute*), 44
Alarm73 (*Alarms attribute*), 44

Alarm74 (*Alarms attribute*), 44
 Alarm75 (*Alarms attribute*), 44
 Alarm76 (*Alarms attribute*), 44
 Alarm77 (*Alarms attribute*), 44
 Alarm78 (*Alarms attribute*), 44
 Alarm79 (*Alarms attribute*), 44
 Alarm8 (*Alarms attribute*), 44
 Alarm8 (*AlarmText attribute*), 40, 63
 Alarm80 (*Alarms attribute*), 44
 Alarm81 (*Alarms attribute*), 44
 Alarm82 (*Alarms attribute*), 44
 Alarm83 (*Alarms attribute*), 44
 Alarm84 (*Alarms attribute*), 44
 Alarm85 (*Alarms attribute*), 44
 Alarm86 (*Alarms attribute*), 44
 Alarm87 (*Alarms attribute*), 44
 Alarm88 (*Alarms attribute*), 44
 Alarm89 (*Alarms attribute*), 44
 Alarm9 (*Alarms attribute*), 44
 Alarm9 (*AlarmText attribute*), 40, 63
 Alarm90 (*Alarms attribute*), 44
 Alarm91 (*Alarms attribute*), 44
 Alarm92 (*Alarms attribute*), 44
 Alarm93 (*Alarms attribute*), 44
 Alarm94 (*Alarms attribute*), 44
 Alarm95 (*Alarms attribute*), 44
 Alarm96 (*Alarms attribute*), 44
 Alarm97 (*Alarms attribute*), 44
 Alarm98 (*Alarms attribute*), 44
 Alarm99 (*Alarms attribute*), 44
 Alarms (*class in hvl_ccb.dev.supercube.constants*), 40
 AlarmText (*class in hvl_ccb.dev.supercube.constants*), 39
 AlarmText (*class in hvl_ccb.dev.supercube2015.constants*), 62
 ALL (*TiePieOscilloscopeAutoResolutionModes attribute*), 82
 all_measurements (*LuminoxMeasurementType attribute*), 170
 all_measurements_types() (*LuminoxMeasurementType class method*), 170
 amplitude (*TiePieGeneratorConfig property*), 78
 ANALOG (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 128
 analog_control (*FuGProbusVDIRegisters property*), 107
 ANY (*LabJack.DeviceType attribute*), 120
 ANY (*LJMCommunicationConfig.ConnectionType attribute*), 16
 ANY (*LJMCommunicationConfig.DeviceType attribute*), 16
 ANY (*TiePieOscilloscopeTriggerKind attribute*), 76
 ARBITRARY (*TiePieGeneratorSignalType attribute*), 79
 AsyncCommunicationProtocol (*class in hvl_ccb.comm.base*), 11
 AsyncCommunicationProtocolConfig (*class in hvl_ccb.comm.base*), 13
 attenuation (*CryLasAttenuator property*), 87
 auto (*EarthingStickOperatingStatus attribute*), 47
 AUTO (*LaudaProRp245eConfig.OperationModeEnum attribute*), 128
 AUTO (*RTO1024.TriggerModes attribute*), 155
 auto_laser_on (*CryLasLaserConfig attribute*), 93
 auto_resolution_mode (*TiePieOscilloscopeConfig property*), 82
 AutoNumberNameEnum (*class in hvl_ccb.utils.enum*), 181

B

B (*SupercubeOpcEndpoint attribute*), 52, 69
 BA (*NewportConfigCommands attribute*), 135
 backlash_compensation (*NewportSMC100PPConfig attribute*), 142
 backup_waveform() (*RTO1024 method*), 155
 bar (*PfeifferTPG.PressureUnits attribute*), 148
 barometric_pressure (*LuminoxMeasurementType attribute*), 170
 base_velocity (*NewportSMC100PPConfig attribute*), 142
 BATH_TEMP (*LaudaProRp245eCommand attribute*), 126
 baudrate (*CryLasAttenuatorSerialCommunicationConfig attribute*), 89
 baudrate (*CryLasLaserSerialCommunicationConfig attribute*), 95
 baudrate (*FuGSerialCommunicationConfig attribute*), 110
 baudrate (*HeinzingerSerialCommunicationConfig attribute*), 116
 baudrate (*LuminoxSerialCommunicationConfig attribute*), 172
 baudrate (*MBW973SerialCommunicationConfig attribute*), 133
 baudrate (*NewportSMC100PPSerialCommunicationConfig attribute*), 146
 baudrate (*PfeifferTPGSerialCommunicationConfig attribute*), 153
 baudrate (*SerialCommunicationConfig attribute*), 24
 BH (*NewportConfigCommands attribute*), 135
 board (*VisaCommunicationConfig attribute*), 31
 BreakdownDetection (*class in hvl_ccb.dev.supercube.constants*), 44
 BreakdownDetection (*class in hvl_ccb.dev.supercube2015.constants*), 63
 bufsize (*TcpCommunicationConfig attribute*), 26
 build_str() (*LaudaProRp245eCommand method*), 127
 bytesize (*CryLasAttenuatorSerialCommunicationConfig attribute*), 89

bytesize (*CryLasLaserSerialCommunicationConfig attribute*), 95
bytesize (*FuGSerialCommunicationConfig attribute*), 110
bytesize (*HeinzingerSerialCommunicationConfig attribute*), 116
bytesize (*LuminoxSerialCommunicationConfig attribute*), 172
bytesize (*MBW973SerialCommunicationConfig attribute*), 133
bytesize (*NewportSMC100PPSerialCommunicationConfig attribute*), 146
bytesize (*PfeifferTPGSerialCommunicationConfig attribute*), 153
Bytesize (*SerialCommunicationConfig attribute*), 24
bytesize (*SerialCommunicationConfig attribute*), 24

C

C (*LabJack.TemperatureUnit attribute*), 120
C (*LabJack.ThermocoupleType attribute*), 120
calibration_factor (*CryLasLaserConfig attribute*), 93
calibration_mode (*FuGProbusVDIRegisters property*), 107
cc_mode (*FuGProbusVDIRegisters property*), 107
cee16 (*GeneralSockets attribute*), 48, 65
channels_enabled (*TiePieOscilloscope property*), 80
check_for_error() (*NewportSMC100PPSerialCommunication method*), 144
check_generic_type() (*in module hvl_ccb.utils.typing*), 182
check_master_slave_config() (*PSI9000 method*), 97
chunk_size (*VisaCommunicationConfig attribute*), 31
clean_amplitude() (*TiePieGeneratorConfig method*), 78
clean_auto_resolution_mode() (*TiePieOscilloscopeConfig static method*), 82
clean_coupling() (*TiePieOscilloscopeChannelConfig static method*), 74
clean_enabled() (*TiePieGeneratorConfig static method*), 78
clean_enabled() (*TiePieOscilloscopeChannelConfig static method*), 74
clean_frequency() (*TiePieGeneratorConfig method*), 78
clean_input_range() (*TiePieOscilloscopeChannelConfig method*), 74
clean_offset() (*TiePieGeneratorConfig method*), 78
clean_pre_sample_ratio() (*TiePieOscilloscopeConfig method*), 82
clean_probe_offset() (*TiePieOscilloscopeChannelConfig method*), 74

clean_record_length() (*TiePieOscilloscopeConfig method*), 82
clean_resolution() (*TiePieOscilloscopeConfig static method*), 82
clean_safe_ground_enabled() (*TiePieOscilloscopeChannelConfig static method*), 74
clean_sample_frequency() (*TiePieOscilloscopeConfig method*), 82
clean_signal_type() (*TiePieGeneratorConfig static method*), 78
clean_trigger_enabled() (*TiePieOscilloscopeChannelConfig static method*), 74
clean_trigger_hysteresis() (*TiePieOscilloscopeChannelConfig method*), 74
clean_trigger_kind() (*TiePieOscilloscopeChannelConfig static method*), 75
clean_trigger_level() (*TiePieOscilloscopeChannelConfig method*), 75
clean_trigger_level_mode() (*TiePieOscilloscopeChannelConfig static method*), 75
clean_trigger_timeout() (*TiePieOscilloscopeConfig method*), 82
clean_values() (*AsyncCommunicationProtocolConfig method*), 13
clean_values() (*CryLasAttenuatorConfig method*), 88
clean_values() (*CryLasLaserConfig method*), 93
clean_values() (*EmptyConfig method*), 85, 184
clean_values() (*FuGConfig method*), 102
clean_values() (*HeinzingerConfig method*), 112
clean_values() (*ILS2TConfig method*), 165
clean_values() (*LaudaProRp245eConfig method*), 128
clean_values() (*LaudaProRp245eTcpCommunicationConfig method*), 130
clean_values() (*LJMCommunicationConfig method*), 17
clean_values() (*LuminoxConfig method*), 169
clean_values() (*MBW973Config method*), 132
clean_values() (*ModbusTcpCommunicationConfig method*), 19
clean_values() (*NewportSMC100PPConfig method*), 142
clean_values() (*OpcUaCommunicationConfig method*), 21
clean_values() (*PfeifferTPGConfig method*), 151
clean_values() (*PSI9000Config method*), 98
clean_values() (*SerialCommunicationConfig method*), 24
clean_values() (*SupercubeConfiguration method*), 37, 60
clean_values() (*TcpCommunicationConfig method*), 26
clean_values() (*TechnixConfig method*), 175

clean_values() (*TelnetCommunicationConfig method*), 29
clean_values() (*TiePieDeviceConfig method*), 72
clean_values() (*VisaCommunicationConfig method*), 31
close (*EarthingStickOperation attribute*), 47
close() (*CommunicationProtocol method*), 14
close() (*LaudaProRp245eTcpCommunication method*), 129
close() (*LJMCommunication method*), 15
close() (*ModbusTcpCommunication method*), 18
close() (*NullCommunicationProtocol method*), 14
close() (*OpcUaCommunication method*), 20
close() (*SerialCommunication method*), 23
close() (*Tcp method*), 26
close() (*TelnetCommunication method*), 28
close() (*VisaCommunication method*), 30
close_shutter() (*CryLasLaser method*), 90
CLOSED (*CryLasLaser.AnswersShutter attribute*), 90
CLOSED (*CryLasLaserShutterStatus attribute*), 96
closed (*DoorStatus attribute*), 45, 63
closed (*EarthingStickStatus attribute*), 47, 65
CMD_EXEC_ERROR (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_CC (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_CONFIGURATION (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_DISABLE (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_HOMING (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_MOVING (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_NOT_REFERENCED (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_PP (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMD_NOT_ALLOWED_READY (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
CMR (*PfeifferTPG.SensorTypes attribute*), 149
CODE_OR_ADDR_INVALID (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
portSMC100PPSerialCommunication.ControllerErrors attribute, 144
collect_measurement_data() (*TiePieOscilloscope method*), 80
com (*SingleCommDevice property*), 86
COM_TIME_OUT (*LaudaProRp245eCommand attribute*), 126
com_time_out (*LaudaProRp245eConfig attribute*), 128
COM_TIMEOUT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 144
command (*LuminoxMeasurementType property*), 170
command() (*FuGProbusIV method*), 105
communication_channel (*TechnixConfig attribute*), 175
CommunicationProtocol (*class in hvl_ccb.comm.base*), 14
config (*ConfigurationMixin property*), 183
CONFIG (*FuGProbusVRegisterGroups attribute*), 108
CONFIG (*NewportSMC100PP.StateMessages attribute*), 136
CONFIG (*NewportStates attribute*), 147
config_cls() (*AsyncCommunicationProtocol static method*), 11
config_cls() (*ConfigurationMixin static method*), 183
config_cls() (*CryLasAttenuator static method*), 87
config_cls() (*CryLasAttenuatorSerialCommunication static method*), 88
config_cls() (*CryLasLaser static method*), 91
config_cls() (*CryLasLaserSerialCommunication static method*), 94
config_cls() (*Device static method*), 84
config_cls() (*FuGProbusIV static method*), 105
config_cls() (*FuGSerialCommunication static method*), 110
config_cls() (*HeinzingerDI static method*), 113
config_cls() (*HeinzingerSerialCommunication static method*), 116
config_cls() (*ILS2T static method*), 163
config_cls() (*ILS2TModbusTcpCommunication static method*), 166
config_cls() (*LaudaProRp245e static method*), 124
config_cls() (*LaudaProRp245eTcpCommunication static method*), 129
config_cls() (*LJMCCommunication static method*), 15
config_cls() (*Luminox static method*), 169
config_cls() (*LuminoxSerialCommunication static method*), 171
config_cls() (*MBW973 static method*), 131
config_cls() (*MBW973SerialCommunication static method*), 18
config_cls() (*NewportSMC100PP static method*), 136
config_cls() (*NewportSMC100PPSerialCommunication*)

static method), 144
config_cls() (*NullCommunicationProtocol static method*), 14
config_cls() (*OpcUaCommunication static method*), 20
config_cls() (*PfeifferTPG static method*), 149
config_cls() (*PfeifferTPGSerialCommunication static method*), 152
config_cls() (*PSI9000 static method*), 97
config_cls() (*PSI9000VisaCommunication static method*), 99
config_cls() (*RTO1024 static method*), 155
config_cls() (*RTO1024VisaCommunication static method*), 161
config_cls() (*SerialCommunication static method*), 23
config_cls() (*Supercube2015Base static method*), 57
config_cls() (*SupercubeAOpcUaCommunication static method*), 53, 71
config_cls() (*SupercubeBase static method*), 33
config_cls() (*SupercubeBOpcUaCommunication static method*), 56
config_cls() (*SupercubeOpcUaCommunication static method*), 38, 61
config_cls() (*SyncCommunicationProtocol static method*), 14
config_cls() (*Tcp static method*), 26
config_cls() (*Technix static method*), 173
config_cls() (*TechnixSerialCommunication static method*), 176
config_cls() (*TechnixTelnetCommunication static method*), 177
config_cls() (*TelnetCommunication static method*), 28
config_cls() (*TiePieOscilloscope static method*), 81
config_cls() (*VisaCommunication static method*), 30
config_cls() (*VisaDevice static method*), 179
config_gen (*TiePieGeneratorMixin attribute*), 78
config_high_pulse () (*LabJack method*), 120
config_i2c (*TiePieI2CHostMixin attribute*), 80
config_osc (*TiePieOscilloscope attribute*), 81
config_osc_channel_dict (*TiePieOscilloscope attribute*), 81
config_status (*FuG property*), 101
configdataclass () (*in module hvl_ccb.configuration*), 184
configuration_save_json () (*ConfigurationMixin method*), 183
ConfigurationMixin (*class in hvl_ccb.configuration*), 183
ConfigurationValueWarning, 183
connection_type (*LJMCommunicationConfig attribute*), 17
CONT_MODE (*LaudaProRp245eCommand attribute*), 126
contact_range () (*GeneralSupport class method*), 48
continue_ramp () (*LaudaProRp245e method*), 124
control_mode (*LaudaProRp245eConfig attribute*), 128
ConvertibleTypes (*in module hvl_ccb.utils.typing*), 182
COOLOFF (*LaudaProRp245eConfig.OperationModeEnum attribute*), 128
COOLON (*LaudaProRp245eConfig.OperationModeEnum attribute*), 128
coupling (*TiePieOscilloscopeChannelConfig property*), 75
CR (*FuGTerminators attribute*), 111
create_serial_port () (*SerialCommunicationConfig method*), 24
create_telnet () (*TelnetCommunicationConfig method*), 29
CRLF (*FuGTerminators attribute*), 111
CryLasAttenuator (*class in hvl_ccb.dev.crylas*), 87
CryLasAttenuatorConfig (*class in hvl_ccb.dev.crylas*), 87
CryLasAttenuatorError, 88
CryLasAttenuatorSerialCommunication (*class in hvl_ccb.dev.crylas*), 88
CryLasAttenuatorSerialCommunicationConfig (*class in hvl_ccb.dev.crylas*), 88
CryLasLaser (*class in hvl_ccb.dev.crylas*), 90
CryLasLaser.AnswersShutter (*class in hvl_ccb.dev.crylas*), 90
CryLasLaser.AnswersStatus (*class in hvl_ccb.dev.crylas*), 90
CryLasLaser.LaserStatus (*class in hvl_ccb.dev.crylas*), 90
CryLasLaser.RepetitionRates (*class in hvl_ccb.dev.crylas*), 90
CryLasLaserConfig (*class in hvl_ccb.dev.crylas*), 93
CryLasLaserError, 93
CryLasLaserNotReadyError, 94
CryLasLaserPoller (*class in hvl_ccb.dev.crylas*), 94
CryLasLaserSerialCommunication (*class in hvl_ccb.dev.crylas*), 94
CryLasLaserSerialCommunicationConfig (*class in hvl_ccb.dev.crylas*), 95
CryLasLaserShutterStatus (*class in hvl_ccb.dev.crylas*), 96
current (*FuG property*), 101
CURRENT (*FuGProbusIVCommands attribute*), 106
CURRENT (*FuGReadbackChannels attribute*), 110
current (*Technix property*), 173
current_lower_limit (*PSI9000Config attribute*), 98
current_monitor (*FuG property*), 101
current_primary (*Power attribute*), 51, 67
current_upper_limit (*PSI9000Config attribute*), 98
CurrentPosition (*New-portSMC100PPConfig.HomeSearch attribute*), 142
cv_mode (*FuGProbusVDIRegisters property*), 107

D

datachange_notification() (*OpcUaSubHandler method*), 22
datachange_notification() (*SupercubeSubscriptionHandler method*), 39, 62
date_of_manufacture (*LuminoxMeasurementType attribute*), 170
DC (*TiePieGeneratorSignalType attribute*), 79
DC_DoubleStage_280kV (*PowerSetup attribute*), 51, 68
DC_SingleStage_140kV (*PowerSetup attribute*), 51, 68
DC_VOLTAGE_TOO_LOW (*NewportSMC100PP.MotorErrors attribute*), 135
DCA (*TiePieOscilloscopeChannelCoupling attribute*), 75
DCV (*TiePieOscilloscopeChannelCoupling attribute*), 75
default_com_cls() (*CryLasAttenuator static method*), 87
default_com_cls() (*CryLasLaser static method*), 91
default_com_cls() (*FuGProbusIV static method*), 105
default_com_cls() (*HeinzingerDI static method*), 113
default_com_cls() (*ILS2T static method*), 163
default_com_cls() (*LabJack static method*), 121
default_com_cls() (*LaudaProRp245e static method*), 125
default_com_cls() (*Luminox static method*), 169
default_com_cls() (*MBW973 static method*), 131
default_com_cls() (*NewportSMC100PP static method*), 136
default_com_cls() (*PfeifferTPG static method*), 149
default_com_cls() (*PSI9000 static method*), 97
default_com_cls() (*RTO1024 static method*), 155
default_com_cls() (*SingleCommDevice static method*), 86
default_com_cls() (*Supercube2015Base static method*), 57
default_com_cls() (*Supercube2015WithFU static method*), 70
default_com_cls() (*SupercubeB static method*), 56
default_com_cls() (*SupercubeBase static method*), 33
default_com_cls() (*SupercubeWithFU static method*), 54
default_com_cls() (*Technix method*), 173
default_com_cls() (*TiePieOscilloscope static method*), 81
default_com_cls() (*VisaDevice static method*), 179
DEFAULT_IO_SCANNING_CONTROL_VALUES (*ILS2T attribute*), 162
default_n_attempts_read_text_nonempty (*AsyncCommunicationProtocolConfig attribute*), 13
default_n_attempts_read_text_nonempty (*FuGSerialCommunicationConfig attribute*), 110
default_n_attempts_read_text_nonempty (*HeinzingerSerialCommunicationConfig attribute*), 116
default_number_of_recordings (*HeinzingerConfig attribute*), 112
Device (*class in hvl_ccb.dev.base*), 84
DEVICE_TYPE (*LaudaProRp245eCommand attribute*), 127
device_type (*LJMCommunicationConfig attribute*), 17
DeviceExistingException, 84
DeviceFailuresException, 84
devices_failed_start (*DeviceSequenceMixin attribute*), 85
devices_failed_start (*ExperimentManager attribute*), 185
devices_failed_stop (*DeviceSequenceMixin attribute*), 85
devices_failed_stop (*ExperimentManager attribute*), 185
DeviceSequenceMixin (*class in hvl_ccb.dev.base*), 85
di (*FuG property*), 101
digital_control (*FuGProbusVDIRegisters property*), 107
DIOChannel (*LabJack attribute*), 119
DISABLE (*NewportStates attribute*), 147
disable() (*ILS2T method*), 163
DISABLE_FROM_JOGGING (*NewportSMC100PP.StateMessages attribute*), 136
DISABLE_FROM_MOVING (*NewportSMC100PP.StateMessages attribute*), 136
DISABLE_FROM_READY (*NewportSMC100PP.StateMessages attribute*), 136
disable_pulses() (*LabJack method*), 121
DISABLED (*TiePieOscilloscopeAutoResolutionModes attribute*), 82
DisableEspStageCheck (*NewportSMC100PPConfig.EspStageConfig attribute*), 142
DISPLACEMENT_OUT_OF_LIMIT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 144
display_message_board() (*SupercubeBase method*), 33
display_status_board() (*SupercubeBase method*), 33
do_ioscanning_write() (*ILS2T method*), 163
Door (*class in hvl_ccb.dev.supercube.constants*), 45
DoorStatus (*class in hvl_ccb.dev.supercube.constants*), 45
DoorStatus (*class in hvl_ccb.dev.supercube2015.constants*), 63

E

E (*LabJack.ThermocoupleType attribute*), 120

E0 (*FuGErrorcodes attribute*), 103
 E1 (*FuGErrorcodes attribute*), 103
 E10 (*FuGErrorcodes attribute*), 103
 E100 (*FuGErrorcodes attribute*), 103
 E106 (*FuGErrorcodes attribute*), 103
 E11 (*FuGErrorcodes attribute*), 103
 E115 (*FuGErrorcodes attribute*), 103
 E12 (*FuGErrorcodes attribute*), 104
 E125 (*FuGErrorcodes attribute*), 104
 E13 (*FuGErrorcodes attribute*), 104
 E135 (*FuGErrorcodes attribute*), 104
 E14 (*FuGErrorcodes attribute*), 104
 E145 (*FuGErrorcodes attribute*), 104
 E15 (*FuGErrorcodes attribute*), 104
 E155 (*FuGErrorcodes attribute*), 104
 E16 (*FuGErrorcodes attribute*), 104
 E165 (*FuGErrorcodes attribute*), 104
 E2 (*FuGErrorcodes attribute*), 104
 E206 (*FuGErrorcodes attribute*), 104
 E306 (*FuGErrorcodes attribute*), 104
 E4 (*FuGErrorcodes attribute*), 104
 E5 (*FuGErrorcodes attribute*), 104
 E504 (*FuGErrorcodes attribute*), 104
 E505 (*FuGErrorcodes attribute*), 104
 E6 (*FuGErrorcodes attribute*), 104
 E666 (*FuGErrorcodes attribute*), 104
 E7 (*FuGErrorcodes attribute*), 104
 E8 (*FuGErrorcodes attribute*), 104
 E9 (*FuGErrorcodes attribute*), 104
EarthingRod (class
 hvl_ccb.dev.supercube.constants), 45
EarthingRodStatus (class
 hvl_ccb.dev.supercube.constants), 45
EarthingStick (class
 hvl_ccb.dev.supercube.constants), 45
EarthingStick (class
 hvl_ccb.dev.supercube2015.constants), 63
EarthingStickMeta (class
 hvl_ccb.dev.supercube.constants), 47
EarthingStickOperatingStatus (class
 hvl_ccb.dev.supercube.constants), 47
EarthingStickOperation (class
 hvl_ccb.dev.supercube.constants), 47
EarthingStickStatus (class
 hvl_ccb.dev.supercube.constants), 47
EarthingStickStatus (class
 hvl_ccb.dev.supercube2015.constants), 65
EEPROM_ACCESS_ERROR (New-
 portSMC100PPSerialCommunication.ControllerErrors
 attribute), 144
EIGHT (*HeinzingerConfig.RecordingsEnum attribute*),
 112
EIGHT_BIT (*TiePieOscilloscopeResolution attribute*), 82
EIGHT_HUNDRED_MILLI_VOLT (*TiePieOscilloscop-
 eRange attribute*), 75
EIGHT_VOLT (*TiePieOscilloscopeRange attribute*), 75
EIGHTBITS (*SerialCommunicationBytesize attribute*), 24
EIGHTY_VOLT (*TiePieOscilloscopeRange attribute*), 75
EmptyConfig (class in *hvl_ccb.configuration*), 183
EmptyConfig (class in *hvl_ccb.dev.base*), 85
enable() (*ILS2T method*), 163
enable_clock() (*LabJack method*), 121
enabled (*TiePieGeneratorConfig property*), 78
enabled (*TiePieOscilloscopeChannelConfig property*),
 75
EnableEspStageCheck (New-
 portSMC100PPConfig.EspStageConfig at-
 tribute), 142
encoding (*AsyncCommunicationProtocolConfig* at-
 tribute), 13
encoding (*NewportSMC100PPSerialCommunicationConfig*
 attribute), 146
encoding_error_handling (*AsyncCommunication-
 ProtocolConfig attribute*), 13
encoding_error_handling (New-
 portSMC100PPSerialCommunicationConfig
 attribute), 146
EndOfRunSwitch (New-
 portSMC100PPConfig.HomeSearch attribute),
 142
EndOfRunSwitch_and_Index (New-
 portSMC100PPConfig.HomeSearch attribute),
 142
in *endpoint_name* (*OpcUaCommunicationConfig* at-
 tribute), 21
in *endpoint_name* (*SupercubeAOpcUaConfiguration* at-
 tribute), 53, 71
in *endpoint_name* (*SupercubeBOpcUaConfiguration* at-
 tribute), 56
in *error* (*DoorStatus attribute*), 45, 63
in *error* (*EarthingStickStatus attribute*), 47, 65
in *ERROR* (*ExperimentStatus attribute*), 186
in *Error* (*SafetyStatus attribute*), 52, 68
in *errorcode* (*FuGError attribute*), 103
in *Errors* (class in *hvl_ccb.dev.supercube.constants*), 48
in *Errors* (class in *hvl_ccb.dev.supercube2015.constants*),
 65
in *ESP_STAGE_NAME_INVALID* (New-
 portSMC100PPSerialCommunication.ControllerErrors
 attribute), 144
in *ETH* (*LaudaProRp245eConfig.ExtControlModeEnum* at-
 tribute), 128
ETHERNET (*LJMCommunicationConfig.ConnectionType*
 attribute), 16
EVEN (*SerialCommunicationParity attribute*), 25
event_notification() (*OpcUaSubHandler method*),
 22

E

- EXECUTE (*FuGProbusIVCommands* attribute), 106
- execute_absolute_position() (*ILS2T* method), 163
- execute_on_x (*FuGProbusVConfigRegisters* property), 107
- execute_relative_step() (*ILS2T* method), 164
- EXECUTEONX (*FuGProbusIVCommands* attribute), 106
- exit_configuration() (*NewportSMC100PP* method), 136
- exit_configuration_wait_sec (*NewportSMC100PPConfig* attribute), 142
- experiment_blocked (*EarthingRodStatus* attribute), 45
- experiment_ready (*EarthingRodStatus* attribute), 45
- ExperimentError, 185
- ExperimentManager (class) in *hvl_ccb.experiment_manager*, 185
- ExperimentStatus (class) in *hvl_ccb.experiment_manager*, 186
- EXPT100 (*LaudaProRp245eConfig*.ExtControlModeEnum attribute), 128
- External (*PowerSetup* attribute), 51, 68
- EXTERNAL_TEMP (*LaudaProRp245eCommand* attribute), 127

F

- F (*LabJack.TemperatureUnit* attribute), 120
- failures (*DeviceFailuresException* attribute), 84
- FALLING (*TiePieOscilloscopeTriggerKind* attribute), 76
- FAST (*ILS2T.Refl6Jog* attribute), 162
- file_copy() (*RTO1024* method), 155
- finish() (*ExperimentManager* method), 185
- FINISHED (*ExperimentStatus* attribute), 186
- FINISHING (*ExperimentStatus* attribute), 186
- FIRMWARE (*FuGRreadbackChannels* attribute), 110
- FIVE_MHZ (*LabJack.ClockFrequency* attribute), 119
- FIVEBITS (*SerialCommunicationBytesize* attribute), 24
- FLT_INFO (*ILS2TRegAddr* attribute), 167
- FLT_MEM_DEL (*ILS2TRegAddr* attribute), 167
- FLT_MEM_RESET (*ILS2TRegAddr* attribute), 167
- FOLLOWING_ERROR (*NewportSMC100PP.MotorErrors* attribute), 135
- FOLLOWRAMP (*FuGRampModes* attribute), 109
- force_trigger() (*TiePieOscilloscope* method), 81
- force_value() (*AsyncCommunicationProtocolConfig* method), 13
- force_value() (*CryLasAttenuatorConfig* method), 88
- force_value() (*CryLasAttenuatorSerialCommunicationConfig* method), 89
- force_value() (*CryLasLaserConfig* method), 93
- force_value() (*CryLasLaserSerialCommunicationConfig* method), 95
- force_value() (*EmptyConfig* method), 86, 184
- force_value() (*FuGConfig* method), 102
- force_value() (*FuGSerialCommunicationConfig* method), 110
- force_value() (*HeinzingerConfig* method), 112
- force_value() (*HeinzingerSerialCommunicationConfig* method), 116
- force_value() (*ILS2TConfig* method), 165
- force_value() (*ILS2TModbusTcpCommunicationConfig* method), 166
- force_value() (*LaudaProRp245eConfig* method), 128
- force_value() (*LaudaProRp245eTcpCommunicationConfig* method), 130
- force_value() (*LJMCommunicationConfig* method), 17
- force_value() (*LuminoxConfig* method), 169
- force_value() (*LuminoxSerialCommunicationConfig* method), 172
- force_value() (*MBW973Config* method), 132
- force_value() (*MBW973SerialCommunicationConfig* method), 133
- force_value() (*ModbusTcpCommunicationConfig* method), 19
- force_value() (*NewportSMC100PPConfig* method), 142
- force_value() (*NewportSMC100PPSerialCommunicationConfig* method), 146
- force_value() (*OpcUaCommunicationConfig* method), 21
- force_value() (*PfeifferTPGConfig* method), 151
- force_value() (*PfeifferTPGSerialCommunicationConfig* method), 153
- force_value() (*PSI9000Config* method), 99
- force_value() (*PSI9000VisaCommunicationConfig* method), 100
- force_value() (*RTO1024Config* method), 160
- force_value() (*RTO1024VisaCommunicationConfig* method), 161
- force_value() (*SerialCommunicationConfig* method), 24
- force_value() (*SupercubeAOpcUaConfiguration* method), 53, 71
- force_value() (*SupercubeBOpcUaConfiguration* method), 56
- force_value() (*SupercubeConfiguration* method), 37, 60
- force_value() (*SupercubeOpcUaCommunicationConfig* method), 38, 61
- force_value() (*TcpCommunicationConfig* method), 27
- force_value() (*TechnixCommunicationConfig* method), 174
- force_value() (*TechnixConfig* method), 175
- force_value() (*TechnixSerialCommunicationConfig* method), 176
- force_value() (*TechnixTelnetCommunicationConfig* method), 177
- force_value() (*TelnetCommunicationConfig* method),

29			
force_value() (<i>TiePieDeviceConfig</i> method), 73			
force_value() (<i>VisaCommunicationConfig</i> method), 31			
force_value() (<i>VisaDeviceConfig</i> method), 180			
FORTY_MHZ (<i>LabJack.ClockFrequency</i> attribute), 119			
FORTY_VOLT (<i>TiePieOscilloscopeRange</i> attribute), 76			
FOUR (<i>HeinzingerConfig.RecordingsEnum</i> attribute), 112			
FOUR_HUNDRED_MILLI_VOLT (<i>TiePieOscilloscopeRange</i> attribute), 76			
FOUR_VOLT (<i>TiePieOscilloscopeRange</i> attribute), 76			
FOURTEEN_BIT (<i>TiePieOscilloscopeResolution</i> attribute), 82			
FREERUN (<i>RTO1024.TriggerModes</i> attribute), 155			
frequency (<i>Power</i> attribute), 51, 67			
frequency (<i>TiePieGeneratorConfig</i> property), 78			
FRM (<i>NewportConfigCommands</i> attribute), 135			
from_json() (<i>ConfigurationMixin</i> class method), 183			
FRS (<i>NewportConfigCommands</i> attribute), 135			
fso_reset() (<i>Supercube2015WithFU</i> method), 70			
fso_reset() (<i>SupercubeWithFU</i> method), 54			
FuG (<i>class in hvl_ccb.dev.fug</i>), 101			
FuGConfig (<i>class in hvl_ccb.dev.fug</i>), 102			
FuDigitalVal (<i>class in hvl_ccb.dev.fug</i>), 103			
FuLError, 103			
FuLErrorcodes (<i>class in hvl_ccb.dev.fug</i>), 103			
FuGMonitorModes (<i>class in hvl_ccb.dev.fug</i>), 104			
FuGPolarities (<i>class in hvl_ccb.dev.fug</i>), 105			
FuGProbusIV (<i>class in hvl_ccb.dev.fug</i>), 105			
FuGProbusIVCommands (<i>class in hvl_ccb.dev.fug</i>), 106			
FuGProbusV (<i>class in hvl_ccb.dev.fug</i>), 106			
FuGProbusVConfigRegisters (<i>class in hvl_ccb.dev.fug</i>), 106			
FuGProbusVDIRegisters (<i>class in hvl_ccb.dev.fug</i>), 107			
FuGProbusVDORegisters (<i>class in hvl_ccb.dev.fug</i>), 108			
FuGProbusVMonitorRegisters (<i>class in hvl_ccb.dev.fug</i>), 108			
FuGProbusVRegisterGroups (<i>class in hvl_ccb.dev.fug</i>), 108			
FuGProbusVSetRegisters (<i>class in hvl_ccb.dev.fug</i>), 109			
FuGRampModes (<i>class in hvl_ccb.dev.fug</i>), 109			
FuGReadbackChannels (<i>class in hvl_ccb.dev.fug</i>), 109			
FuGSerialCommunication (<i>class in hvl_ccb.dev.fug</i>), 110			
FuGSerialCommunicationConfig (<i>class in hvl_ccb.dev.fug</i>), 110			
FuGTerminators (<i>class in hvl_ccb.dev.fug</i>), 111			
G			
GeneralSockets (<i>class in hvl_ccb.dev.supercube.constants</i>), 48			
GeneralSockets (<i>class in hvl_ccb.dev.supercube2015.constants</i>), 65			
GeneralSupport (<i>class in hvl_ccb.dev.supercube.constants</i>), 48			
GeneralSupport (<i>class in hvl_ccb.dev.supercube2015.constants</i>), 65			
GeneralSupportMeta (<i>class in hvl_ccb.dev.supercube.constants</i>), 49			
GENERATOR (<i>TiePieDeviceType</i> attribute), 73			
generator_start() (<i>TiePieGeneratorMixin</i> method), 78			
generator_stop() (<i>TiePieGeneratorMixin</i> method), 79			
get() (<i>AlarmText</i> class method), 40, 63			
get() (<i>MeasurementsDividerRatio</i> class method), 67			
get() (<i>MeasurementsScaledInput</i> class method), 67			
get_acceleration() (<i>NewportSMC100PP</i> method), 137			
get_acquire_length() (<i>RTO1024</i> method), 155			
get_ain() (<i>LabJack</i> method), 121			
get_bath_temp() (<i>LaudaProRp245e</i> method), 125			
get_by_p_id() (<i>LabJack.DeviceType</i> class method), 120			
get_by_p_id() (<i>LJMCommunicationConfig.DeviceType</i> class method), 17			
get_cal_current_source() (<i>LabJack</i> method), 121			
get_cee16_socket() (<i>Supercube2015Base</i> method), 57			
get_cee16_socket() (<i>SupercubeBase</i> method), 33			
get_channel_offset() (<i>RTO1024</i> method), 156			
get_channel_position() (<i>RTO1024</i> method), 156			
get_channel_range() (<i>RTO1024</i> method), 156			
get_channel_scale() (<i>RTO1024</i> method), 156			
get_channel_state() (<i>RTO1024</i> method), 156			
get_clock() (<i>LabJack</i> method), 121			
get_controller_information() (<i>NewportSMC100PP</i> method), 137			
get_current() (<i>HeinzingerDI</i> method), 113			
get_dc_volt() (<i>ILS2T</i> method), 164			
get_device() (<i>DeviceSequenceMixin</i> method), 85			
get_device_by_serial_number() (<i>in module hvl_ccb.dev.tiepie.base</i>), 73			
get_device_type() (<i>LaudaProRp245e</i> method), 125			
get_devices() (<i>DeviceSequenceMixin</i> method), 85			
get_digital_input() (<i>LabJack</i> method), 121			
get_door_status() (<i>Supercube2015Base</i> method), 57			
get_door_status() (<i>SupercubeBase</i> method), 33			
get_earthing_manual() (<i>Supercube2015Base</i> method), 57			
get_earthing_rod_status() (<i>SupercubeBase</i> method), 34			
get_earthing_status() (<i>Supercube2015Base</i> method), 58			
get_earthing_stick_manual() (<i>SupercubeBase</i> method), 34			

get_earthing_stick_operating_status() (*SupercubeBase method*), 34
 get_earthing_stick_status() (*SupercubeBase method*), 34
 get_error_code() (*ILS2T method*), 164
 get_error_queue() (*VisaDevice method*), 179
 get_frequency() (*Supercube2015WithFU method*), 70
 get_frequency() (*SupercubeWithFU method*), 54
 get_fso_active() (*Supercube2015WithFU method*), 70
 get_fso_active() (*SupercubeWithFU method*), 54
 get_full_scale_mbar() (*PfeifferTPG method*), 149
 get_full_scale_unitless() (*PfeifferTPG method*), 149
 get_identification() (*VisaDevice method*), 179
 get_interface_version() (*HeinzingerDI method*), 113
 get_max_voltage() (*Supercube2015WithFU method*), 70
 get_max_voltage() (*SupercubeWithFU method*), 54
 get_measurement_ratio() (*Supercube2015Base method*), 58
 get_measurement_ratio() (*SupercubeBase method*), 34
 get_measurement_voltage() (*Supercube2015Base method*), 58
 get_measurement_voltage() (*SupercubeBase method*), 34
 get_motor_configuration() (*NewportSMC100PP method*), 137
 get_move_duration() (*NewportSMC100PP method*), 137
 get_negative_software_limit() (*NewportSMC100PP method*), 138
 get_number_of_recordings() (*HeinzingerDI method*), 113
 get_output() (*PSI9000 method*), 97
 get_position() (*ILS2T method*), 164
 get_position() (*NewportSMC100PP method*), 138
 get_positive_software_limit() (*NewportSMC100PP method*), 138
 get_power_setup() (*Supercube2015WithFU method*), 70
 get_power_setup() (*SupercubeWithFU method*), 54
 get_primary_current() (*Supercube2015WithFU method*), 71
 get_primary_current() (*SupercubeWithFU method*), 54
 get_primary_voltage() (*Supercube2015WithFU method*), 71
 get_primary_voltage() (*SupercubeWithFU method*), 54
 get_product_id() (*LabJack method*), 121
 get_product_name() (*LabJack method*), 121
 get_product_type() (*LabJack method*), 122
 get_pulse_energy_and_rate() (*CryLasLaser method*), 91
 get_reference_point() (*RTO1024 method*), 156
 get_register() (*FuGProbusV method*), 106
 get_repetitions() (*RTO1024 method*), 156
 get_sbus_rh() (*LabJack method*), 122
 get_sbus_temp() (*LabJack method*), 122
 get_serial_number() (*HeinzingerDI method*), 113
 get_serial_number() (*LabJack method*), 122
 get_state() (*NewportSMC100PP method*), 138
 get_status() (*ILS2T method*), 164
 get_status() (*Supercube2015Base method*), 58
 get_status() (*SupercubeBase method*), 34
 get_status_byte() (*Technix method*), 173
 get_support_input() (*Supercube2015Base method*), 58
 get_support_input() (*SupercubeBase method*), 34
 get_support_output() (*Supercube2015Base method*), 58
 get_support_output() (*SupercubeBase method*), 35
 get_system_lock() (*PSI9000 method*), 97
 get_t13_socket() (*Supercube2015Base method*), 58
 get_t13_socket() (*SupercubeBase method*), 35
 get_target_voltage() (*Supercube2015WithFU method*), 71
 get_target_voltage() (*SupercubeWithFU method*), 55
 get_temperature() (*ILS2T method*), 164
 get_timestamps() (*RTO1024 method*), 156
 get_ui_lower_limits() (*PSI9000 method*), 97
 get_uip_upper_limits() (*PSI9000 method*), 97
 get_voltage() (*HeinzingerDI method*), 114
 get_voltage_current_setpoint() (*PSI9000 method*), 97
 go_home() (*NewportSMC100PP method*), 138
 go_to_configuration() (*NewportSMC100PP method*), 139
 GreenNotReady (*SafetyStatus attribute*), 52, 69
 GreenReady (*SafetyStatus attribute*), 52, 69

H

HARDWARE (*CryLasLaser.RepetitionRates attribute*), 90
 has_safe_ground (*TiePieOscilloscopeChannelConfig property*), 75
 HEAD (*CryLasLaser.AnswersStatus attribute*), 90
 HeinzingerConfig (*class in hvl_ccb.dev.heinzinger*), 112
 HeinzingerConfig.RecordingsEnum (*class in hvl_ccb.dev.heinzinger*), 112
 HeinzingerDI (*class in hvl_ccb.dev.heinzinger*), 113
 HeinzingerDI.OutputStatus (*class in hvl_ccb.dev.heinzinger*), 113
 HeinzingerPNC (*class in hvl_ccb.dev.heinzinger*), 115

HeinzingerPNC.UnitCurrent	(class <i>hvl_ccb.dev.heinzinger</i>), 115	<i>in</i>	module, 11
HeinzingerPNC.UnitVoltage	(class <i>hvl_ccb.dev.heinzinger</i>), 115	<i>in</i>	<i>hvl_ccb.comm.labjack_ljm</i> module, 15
HeinzingerPNCDeviceNotRecognizedException,	115		<i>hvl_ccb.comm.modbus_tcp</i> module, 18
HeinzingerPNCError	116		<i>hvl_ccb.comm.opc</i> module, 20
HeinzingerPNCMaxCurrentExceededException	116		<i>hvl_ccb.comm.serial</i> module, 23
HeinzingerPNCMaxVoltageExceededException	116		<i>hvl_ccb.comm.tcp</i> module, 26
HeinzingerSerialCommunication	(class <i>hvl_ccb.dev.heinzinger</i>), 116	<i>in</i>	<i>hvl_ccb.comm.telnet</i> module, 28
HeinzingerSerialCommunicationConfig	(class <i>hvl_ccb.dev.heinzinger</i>), 116	<i>in</i>	<i>hvl_ccb.comm.visa</i> module, 30
HIGH (<i>LabJack.DIOStatus</i> attribute)	119		<i>hvl_ccb.configuration</i> module, 183
high_resolution (<i>FuGProbusVSetRegisters</i> property),	109		<i>hvl_ccb.dev</i> module, 181
home_search_polling_interval	(<i>NewportSMC100PPConfig</i> attribute), 143		<i>hvl_ccb.dev.base</i> module, 84
home_search_timeout (<i>NewportSMC100PPConfig</i> attribute)	143		<i>hvl_ccb.dev.crylas</i> module, 87
home_search_type (<i>NewportSMC100PPConfig</i> attribute)	143		<i>hvl_ccb.dev.ea_psi9000</i> module, 96
home_search_velocity (<i>NewportSMC100PPConfig</i> attribute)	143		<i>hvl_ccb.dev.fug</i> module, 100
HOME_STARTED (<i>NewportSMC100PPSerialCommunication</i> . <i>Cont</i> attribute), 144			<i>hvl_ccb.dev.heinzinger</i> module, 112
HomeSwitch	(<i>NewportSMC100PPConfig.HomeSearch</i> attribute), 142		<i>hvl_ccb.dev.labjack</i> module, 118
HomeSwitch_and_Index	(<i>NewportSMC100PPConfig.HomeSearch</i> attribute), 142		<i>hvl_ccb.dev.lauda</i> module, 124
HOMING (<i>NewportStates</i> attribute)	147		<i>hvl_ccb.dev.mbw973</i> module, 131
HOMING_FROM_RS232	(<i>NewportSMC100PP.StateMessages</i> attribute), 136		<i>hvl_ccb.dev.newport</i> module, 134
HOMING_FROM_SMC	(<i>NewportSMC100PP.StateMessages</i> attribute), 136		<i>hvl_ccb.dev.pfeiffer_tpg</i> module, 148
HOMING_TIMEOUT (<i>NewportSMC100PP.MotorErrors</i> attribute)	135		<i>hvl_ccb.dev.rs_rto1024</i> module, 154
horn (<i>Safety</i> attribute)	68		<i>hvl_ccb.dev.se_ilst2t</i> module, 162
horn() (<i>Supercube2015Base</i> method)	59		<i>hvl_ccb.dev.sst_luminox</i> module, 168
host (<i>ModbusTcpCommunicationConfig</i> attribute)	19		<i>hvl_ccb.dev.supercube</i> module, 57
host (<i>OpcUaCommunicationConfig</i> attribute)	21		<i>hvl_ccb.dev.supercube.base</i> module, 33
host (<i>TcpCommunicationConfig</i> attribute)	27		<i>hvl_ccb.dev.supercube.constants</i> module, 39
host (<i>TelnetCommunicationConfig</i> attribute)	29		<i>hvl_ccb.dev.supercube.typ_a</i> module, 53
host (<i>VisaCommunicationConfig</i> attribute)	32		<i>hvl_ccb.dev.supercube.typ_b</i>
hPascal (<i>PfeifferTPG.PressureUnits</i> attribute)	148		
HT (<i>NewportConfigCommands</i> attribute)	135		
hv (<i>Technix</i> property)	173		
hvl_ccb			
<i>module</i> , 186			
hvl_ccb.comm			
<i>module</i> , 33			
hvl_ccb.comm.base			

module, 55
hvl_ccb.dev.supercube2015
 module, 72
hvl_ccb.dev.supercube2015.base
 module, 57
hvl_ccb.dev.supercube2015.constants
 module, 62
hvl_ccb.dev.supercube2015.typ_a
 module, 70
hvl_ccb.dev.technix
 module, 173
hvl_ccb.dev.tiepie
 module, 83
hvl_ccb.dev.tiepie.base
 module, 72
hvl_ccb.dev.tiepie.channel
 module, 74
hvl_ccb.dev.tiepie.device
 module, 77
hvl_ccb.dev.tiepie.generator
 module, 78
hvl_ccb.dev.tiepie.i2c
 module, 79
hvl_ccb.dev.tiepie.oscilloscope
 module, 80
hvl_ccb.dev.tiepie.utils
 module, 83
hvl_ccb.dev.utils
 module, 177
hvl_ccb.dev.visa
 module, 179
hvl_ccb.experiment_manager
 module, 185
hvl_ccb.utils
 module, 183
hvl_ccb.utils.enum
 module, 181
hvl_ccb.utils.typing
 module, 182
hvl_ccb.utils.validation
 module, 182
hysteresis_compensation (New)
portSMC100PPConfig attribute, 143

|

I2C (*TiePieDeviceType attribute*), 73
ID (*FuGProbusIVCommands attribute*), 106
Identification_error (*PfeifferTPG.SensorStatus attribute*), 148
identifier (*LJMCommunicationConfig attribute*), 17
identify_device() (*FuG method*), 101
identify_device() (*HeinzingerPNC method*), 115
identify_sensors() (*PfeifferTPG method*), 149
IKR (*PfeifferTPG.SensorTypes attribute*), 149

IKR11 (*PfeifferTPG.SensorTypes attribute*), 149
IKR9 (*PfeifferTPG.SensorTypes attribute*), 149
ILS2T (*class in hvl_ccb.dev.se_il2t*), 162
ILS2T.ActionsPtp (*class in hvl_ccb.dev.se_il2t*), 162
ILS2T.Mode (*class in hvl_ccb.dev.se_il2t*), 162
ILS2T.Ref16Jog (*class in hvl_ccb.dev.se_il2t*), 162
ILS2T.State (*class in hvl_ccb.dev.se_il2t*), 163
ILS2TConfig (*class in hvl_ccb.dev.se_il2t*), 165
ILS2TException, 166
ILS2TModbusTcpCommunication (*class in hvl_ccb.dev.se_il2t*), 166
ILS2TModbusTcpCommunicationConfig (*class in hvl_ccb.dev.se_il2t*), 166
ILS2TRegAddr (*class in hvl_ccb.dev.se_il2t*), 167
ILS2TRegDatatype (*class in hvl_ccb.dev.se_il2t*), 167
IMMEDIATELY (*FuGRampModes attribute*), 109
IMR (*PfeifferTPG.SensorTypes attribute*), 149
in_1_1 (*GeneralSupport attribute*), 48, 65
in_1_2 (*GeneralSupport attribute*), 48, 65
in_2_1 (*GeneralSupport attribute*), 48, 66
in_2_2 (*GeneralSupport attribute*), 48, 66
in_3_1 (*GeneralSupport attribute*), 48, 66
in_3_2 (*GeneralSupport attribute*), 48, 66
in_4_1 (*GeneralSupport attribute*), 48, 66
in_4_2 (*GeneralSupport attribute*), 48, 66
in_5_1 (*GeneralSupport attribute*), 48, 66
in_5_2 (*GeneralSupport attribute*), 48, 66
in_6_1 (*GeneralSupport attribute*), 48, 66
in_6_2 (*GeneralSupport attribute*), 49, 66
INACTIVE (*CryLasLaser.AnswersStatus attribute*), 90
inactive (*DoorStatus attribute*), 45, 63
inactive (*EarthingStickStatus attribute*), 47, 65
inhibit (*Technix property*), 173
init_attenuation (*CryLasAttenuatorConfig attribute*), 88
init_monitored_nodes() (*OpcUaCommunication method*), 20
init_shutter_status (*CryLasLaserConfig attribute*), 93
initialize() (*NewportSMC100PP method*), 139
INITIALIZED (*ExperimentStatus attribute*), 186
INITIALIZING (*ExperimentStatus attribute*), 186
Initializing (*SafetyStatus attribute*), 52, 69
INPUT (*FuGProbusVRegisterGroups attribute*), 108
input() (*GeneralSupport class method*), 49, 66
input_1 (*MeasurementsDividerRatio attribute*), 49, 67
input_1 (*MeasurementsScaledInput attribute*), 50, 67
input_2 (*MeasurementsDividerRatio attribute*), 50
input_2 (*MeasurementsScaledInput attribute*), 50, 67
input_3 (*MeasurementsDividerRatio attribute*), 50
input_3 (*MeasurementsScaledInput attribute*), 50, 67
input_4 (*MeasurementsDividerRatio attribute*), 50
input_4 (*MeasurementsScaledInput attribute*), 50, 67

input_range (*TiePieOscilloscopeChannelConfig* property), 75
 INT32 (*ILS2TRegDatatype* attribute), 168
 interface_type (*PSI9000VisaCommunicationConfig* attribute), 100
 interface_type (*RTO1024VisaCommunicationConfig* attribute), 161
 interface_type (*VisaCommunicationConfig* attribute), 32
 internal (*LabJack.CjcType* attribute), 119
 INTERNAL (*LaudaProRp245eConfig.ExtControlModeEnum* attribute), 128
 Internal (*PowerSetup* attribute), 51, 68
 InvalidSupercubeStatusError, 57
 IO_SCANNING (*ILS2TRegAddr* attribute), 167
 IoScanningModeValueError, 168
 is_configdataclass (*AsyncCommunicationProtocolConfig* attribute), 13
 is_configdataclass (*CryLasAttenuatorConfig* attribute), 88
 is_configdataclass (*CryLasLaserConfig* attribute), 93
 is_configdataclass (*EmptyConfig* attribute), 86, 184
 is_configdataclass (*FuGConfig* attribute), 102
 is_configdataclass (*HeinzingerConfig* attribute), 112
 is_configdataclass (*ILS2TConfig* attribute), 166
 is_configdataclass (*LaudaProRp245eConfig* attribute), 128
 is_configdataclass (*LJMCommunicationConfig* attribute), 17
 is_configdataclass (*LuminoxConfig* attribute), 170
 is_configdataclass (*MBW973Config* attribute), 132
 is_configdataclass (*ModbusTcpCommunicationConfig* attribute), 19
 is_configdataclass (*NewportSMC100PPConfig* attribute), 143
 is_configdataclass (*OpcUaCommunicationConfig* attribute), 21
 is_configdataclass (*PfeifferTPGConfig* attribute), 151
 is_configdataclass (*SupercubeConfiguration* attribute), 37, 60
 is_configdataclass (*TcpCommunicationConfig* attribute), 27
 is_configdataclass (*TechnixConfig* attribute), 175
 is_configdataclass (*TiePieDeviceConfig* attribute), 73
 is_configdataclass (*VisaCommunicationConfig* attribute), 32
 is_data_ready_polling_interval_sec (*TiePieDeviceConfig* attribute), 73
 is_done() (*MBW973* method), 131
 is_error() (*ExperimentManager* method), 185
 is_finished() (*ExperimentManager* method), 186
 is_generic_type_hint() (in *hvl_ccb.utils.typing*), 182
 is_in_range() (*ILS2TRegDatatype* method), 168
 is_inactive (*CryLasLaser.LaserStatus* property), 90
 is_measurement_data_ready() (*TiePieOscilloscope* method), 81
 is_open (*LJMCommunication* property), 15
 is_open (*OpcUaCommunication* property), 20
 is_open (*SerialCommunication* property), 23
 is_open (*TelnetCommunication* property), 28
 is_polling() (*Poller* method), 178
 is_ready (*CryLasLaser.LaserStatus* property), 90
 is_running() (*ExperimentManager* method), 186
 is_triggered() (*TiePieOscilloscope* method), 81
 is_valid_scale_range_reversed_str() (*PfeifferTPGConfig.Model* method), 151

J

J (*LabJack.ThermocoupleType* attribute), 120
 jerk_time (*NewportSMC100PPConfig* attribute), 143
 JOG (*ILS2T.Mode* attribute), 162
 jog_run() (*ILS2T* method), 164
 jog_stop() (*ILS2T* method), 164
 JOGGING (*NewportStates* attribute), 147
 JOGGING_FROM_DISABLE (*NewportSMC100PP.StateMessages* attribute), 136
 JOGGING_FROM_READY (*NewportSMC100PP.StateMessages* attribute), 136
 JOGN_FAST (*ILS2TRegAddr* attribute), 167
 JOGN_SLOW (*ILS2TRegAddr* attribute), 167
 JR (*NewportConfigCommands* attribute), 135

K

K (*LabJack.TemperatureUnit* attribute), 120
 K (*LabJack.ThermocoupleType* attribute), 120
 keys() (*AsyncCommunicationProtocolConfig* class method), 13
 keys() (*CryLasAttenuatorConfig* class method), 88
 keys() (*CryLasAttenuatorSerialCommunicationConfig* class method), 89
 keys() (*CryLasLaserConfig* class method), 93
 keys() (*CryLasLaserSerialCommunicationConfig* class method), 95
 keys() (*EmptyConfig* class method), 86, 184
 keys() (*FuGConfig* class method), 102
 keys() (*FuGSerialCommunicationConfig* class method), 111
 keys() (*HeinzingerConfig* class method), 112
 keys() (*HeinzingerSerialCommunicationConfig* class method), 117
 keys() (*ILS2TConfig* class method), 166

<code>keys()</code> (<i>ILS2TModbusTcpCommunicationConfig class method</i>), 167	<code>LabJack.CalMicroAmpere</code> (class in <i>hvl_ccb.dev.labjack</i>), 119
<code>keys()</code> (<i>LaudaProRp245eConfig class method</i>), 129	<code>LabJack.CjcType</code> (class in <i>hvl_ccb.dev.labjack</i>), 119
<code>keys()</code> (<i>LaudaProRp245eTcpCommunicationConfig class method</i>), 130	<code>LabJack.ClockFrequency</code> (class in <i>hvl_ccb.dev.labjack</i>), 119
<code>keys()</code> (<i>LJMCommunicationConfig class method</i>), 17	<code>LabJack.DeviceType</code> (class in <i>hvl_ccb.dev.labjack</i>), 119
<code>keys()</code> (<i>LuminoxConfig class method</i>), 170	<code>LabJack.DIOStatus</code> (class in <i>hvl_ccb.dev.labjack</i>), 119
<code>keys()</code> (<i>LuminoxSerialCommunicationConfig class method</i>), 172	<code>LabJack.TemperatureUnit</code> (class in <i>hvl_ccb.dev.labjack</i>), 120
<code>keys()</code> (<i>MBW973Config class method</i>), 132	<code>LabJack.ThermocoupleType</code> (class in <i>hvl_ccb.dev.labjack</i>), 120
<code>keys()</code> (<i>MBW973SerialCommunicationConfig class method</i>), 134	<code>LabJackError</code> , 124
<code>keys()</code> (<i>ModbusTcpCommunicationConfig class method</i>), 19	<code>LabJackIdentifierDIOError</code> , 124
<code>keys()</code> (<i>NewportSMC100PPConfig class method</i>), 143	<code>laser_off()</code> (<i>CryLasLaser method</i>), 91
<code>keys()</code> (<i>NewportSMC100PPSerialCommunicationConfig class method</i>), 146	<code>laser_on()</code> (<i>CryLasLaser method</i>), 91
<code>keys()</code> (<i>OpcUaCommunicationConfig class method</i>), 21	<code>LaudaProRp245e</code> (class in <i>hvl_ccb.dev.lauda</i>), 124
<code>keys()</code> (<i>PfeifferTPGConfig class method</i>), 151	<code>LaudaProRp245eCommand</code> (class in <i>hvl_ccb.dev.lauda</i>), 126
<code>keys()</code> (<i>PfeifferTPGSerialCommunicationConfig class method</i>), 153	<code>LaudaProRp245eCommandError</code> , 127
<code>keys()</code> (<i>PSI9000Config class method</i>), 99	<code>LaudaProRp245eConfig</code> (class in <i>hvl_ccb.dev.lauda</i>), 128
<code>keys()</code> (<i>PSI9000VisaCommunicationConfig class method</i>), 100	<code>LaudaProRp245eConfig.ExtControlModeEnum</code> (class in <i>hvl_ccb.dev.lauda</i>), 128
<code>keys()</code> (<i>RTO1024Config class method</i>), 160	<code>LaudaProRp245eConfig.OperationModeEnum</code> (class in <i>hvl_ccb.dev.lauda</i>), 128
<code>keys()</code> (<i>RTO1024VisaCommunicationConfig class method</i>), 161	<code>LaudaProRp245eTcpCommunication</code> (class in <i>hvl_ccb.dev.lauda</i>), 129
<code>keys()</code> (<i>SerialCommunicationConfig class method</i>), 25	<code>LaudaProRp245eTcpCommunicationConfig</code> (class in <i>hvl_ccb.dev.lauda</i>), 130
<code>keys()</code> (<i>SupercubeAOpcUaConfiguration class method</i>), 53, 71	<code>LF</code> (<i>FuGTerminators attribute</i>), 111
<code>keys()</code> (<i>SupercubeBOpcUaConfiguration class method</i>), 56	<code>LFCR</code> (<i>FuGTerminators attribute</i>), 111
<code>keys()</code> (<i>SupercubeConfiguration class method</i>), 37, 60	<code>line_1</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>SupercubeOpcUaCommunicationConfig class method</i>), 38, 61	<code>line_10</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TcpCommunicationConfig class method</i>), 27	<code>line_11</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TechnixCommunicationConfig class method</i>), 174	<code>line_12</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TechnixConfig class method</i>), 175	<code>line_13</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TechnixSerialCommunicationConfig class method</i>), 176	<code>line_14</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TechnixTelnetCommunicationConfig class method</i>), 177	<code>line_15</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TelnetCommunicationConfig class method</i>), 29	<code>line_2</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>TiePieDeviceConfig class method</i>), 73	<code>line_3</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>VisaCommunicationConfig class method</i>), 32	<code>line_4</code> (<i>MessageBoard attribute</i>), 50
<code>keys()</code> (<i>VisaDeviceConfig class method</i>), 180	<code>line_5</code> (<i>MessageBoard attribute</i>), 50
<code>kV</code> (<i>HeinzingerPNC.UnitVoltage attribute</i>), 115	<code>line_6</code> (<i>MessageBoard attribute</i>), 50
L	<code>line_7</code> (<i>MessageBoard attribute</i>), 50
<code>LabJack</code> (class in <i>hvl_ccb.dev.labjack</i>), 118	<code>line_8</code> (<i>MessageBoard attribute</i>), 50
<code>LabJack.AInRange</code> (class in <i>hvl_ccb.dev.labjack</i>), 118	<code>line_9</code> (<i>MessageBoard attribute</i>), 50
<code>LabJack.BitLimit</code> (class in <i>hvl_ccb.dev.labjack</i>), 119	<code>list_devices()</code> (<i>TiePieOscilloscope static method</i>), 81
	<code>list_directory()</code> (<i>RTO1024 method</i>), 156
	<code>live</code> (<i>OpcControl attribute</i>), 50
	<code>LJMCommunication</code> (class in <i>hvl_ccb.comm.labjack_ljm</i>), 15
	<code>LJMCommunicationConfig</code> (class in <i>hvl_ccb.comm.labjack_ljm</i>), 16

LJMCommunicationConfig.ConnectionType (class in *hvl_ccb.comm.labjack_ljm*), 16
 LJMCommunicationConfig.DeviceType (class in *hvl_ccb.comm.labjack_ljm*), 16
 LJMCommunicationError, 17
 lm34 (*LabJack.CjcType* attribute), 119
 load_configuration() (*RTO1024* method), 157
 local_display() (*RTO1024* method), 157
 locked (*DoorStatus* attribute), 45, 63
 LOW (*LabJack.DIOStatus* attribute), 119
 LOWER_TEMP (*LaudaProRp245eCommand* attribute), 127
 lower_temp (*LaudaProRp245eConfig* attribute), 129
 Luminox (class in *hvl_ccb.dev.sst_luminox*), 168
 LuminoxConfig (class in *hvl_ccb.dev.sst_luminox*), 169
 LuminoxMeasurementType (class in *hvl_ccb.dev.sst_luminox*), 170
 LuminoxMeasurementTypeDict (in module *hvl_ccb.dev.sst_luminox*), 171
 LuminoxMeasurementTypeError, 171
 LuminoxMeasurementTypeValue (in module *hvl_ccb.dev.sst_luminox*), 171
 LuminoxOutputMode (class in *hvl_ccb.dev.sst_luminox*), 171
 LuminoxOutputModeError, 171
 LuminoxSerialCommunication (class in *hvl_ccb.dev.sst_luminox*), 171
 LuminoxSerialCommunicationConfig (class in *hvl_ccb.dev.sst_luminox*), 171

M

mA (*HeinzingerPNC.UnitCurrent* attribute), 115
 manual (*EarthingStickOperatingStatus* attribute), 47
 manual() (*EarthingStick* class method), 46, 64
 manual_1 (*EarthingStick* attribute), 46, 64
 manual_2 (*EarthingStick* attribute), 46, 64
 manual_3 (*EarthingStick* attribute), 46, 64
 manual_4 (*EarthingStick* attribute), 46, 64
 manual_5 (*EarthingStick* attribute), 46, 64
 manual_6 (*EarthingStick* attribute), 46, 64
 manuals() (*EarthingStick* class method), 46
 MARK (*SerialCommunicationParity* attribute), 25
 max_current (*FuG* property), 101
 max_current (*HeinzingerPNC* property), 115
 max_current (*Technix* property), 173
 max_current (*TechnixConfig* attribute), 175
 max_current_hardware (*FuG* property), 101
 max_current_hardware (*HeinzingerPNC* property), 115
 max_pr_number (*LaudaProRp245eConfig* attribute), 129
 max_pump_level (*LaudaProRp245eConfig* attribute), 129
 max_timeout_retry_nr (*OpcUaCommunicationConfig* attribute), 21
 max_voltage (*FuG* property), 101

max_voltage (*HeinzingerPNC* property), 115
 max_voltage (*Technix* property), 173
 max_voltage (*TechnixConfig* attribute), 175
 max_voltage_hardware (*FuG* property), 101
 max_voltage_hardware (*HeinzingerPNC* property), 115
 MAXIMUM (*LabJack.ClockFrequency* attribute), 119
 mbar (*PfeifferTPG.PressureUnits* attribute), 148
 MBW973 (class in *hvl_ccb.dev.mbw973*), 131
 MBW973Config (class in *hvl_ccb.dev.mbw973*), 132
 MBW973ControlRunningException, 133
 MBW973Error, 133
 MBW973PumpRunningException, 133
 MBW973SerialCommunication (class in *hvl_ccb.dev.mbw973*), 133
 MBW973SerialCommunicationConfig (class in *hvl_ccb.dev.mbw973*), 133
 measure() (*PfeifferTPG* method), 150
 measure_all() (*PfeifferTPG* method), 150
 measure_current() (*HeinzingerDI* method), 114
 measure_voltage() (*HeinzingerDI* method), 114
 measure_voltage_current() (*PSI9000* method), 97
 MeasurementsDividerRatio (class in *hvl_ccb.dev.supercube.constants*), 49
 MeasurementsDividerRatio (class in *hvl_ccb.dev.supercube2015.constants*), 66
 MeasurementsScaledInput (class in *hvl_ccb.dev.supercube.constants*), 50
 MeasurementsScaledInput (class in *hvl_ccb.dev.supercube2015.constants*), 67
 message (*Errors* attribute), 48
 MessageBoard (class in *hvl_ccb.dev.supercube.constants*), 50
 micro_step_per_full_step_factor (New-portSMC100PPConfig attribute), 143
 Micron (*PfeifferTPG.PressureUnits* attribute), 148
 MINIMUM (*LabJack.ClockFrequency* attribute), 119
 ModbusTcpCommunication (class in *hvl_ccb.comm.modbus_tcp*), 18
 ModbusTcpCommunicationConfig (class in *hvl_ccb.comm.modbus_tcp*), 19
 ModbusTcpConnectionFailedException, 19
 model (*PfeifferTPGConfig* attribute), 151
 module
 hvl_ccb, 186
 hvl_ccb.comm, 33
 hvl_ccb.comm.base, 11
 hvl_ccb.comm.labjack_ljm, 15
 hvl_ccb.comm.modbus_tcp, 18
 hvl_ccb.comm.opc, 20
 hvl_ccb.comm.serial, 23
 hvl_ccb.comm.tcp, 26
 hvl_ccb.comm.telnet, 28
 hvl_ccb.comm.visa, 30

hvl_ccb.configuration, 183
 hvl_ccb.dev, 181
 hvl_ccb.dev.base, 84
 hvl_ccb.dev.crylas, 87
 hvl_ccb.dev.ea_psi9000, 96
 hvl_ccb.dev.fug, 100
 hvl_ccb.dev.heinzinger, 112
 hvl_ccb.dev.labjack, 118
 hvl_ccb.dev.lauda, 124
 hvl_ccb.dev.mbw973, 131
 hvl_ccb.dev.newport, 134
 hvl_ccb.dev.pfeiffer_tpg, 148
 hvl_ccb.dev.rs_rto1024, 154
 hvl_ccb.dev.se_ilst2t, 162
 hvl_ccb.dev.sst_luminox, 168
 hvl_ccb.dev.supercube, 57
 hvl_ccb.dev.supercube.base, 33
 hvl_ccb.dev.supercube.constants, 39
 hvl_ccb.dev.supercube.typ_a, 53
 hvl_ccb.dev.supercube.typ_b, 55
 hvl_ccb.dev.supercube2015, 72
 hvl_ccb.dev.supercube2015.base, 57
 hvl_ccb.dev.supercube2015.constants, 62
 hvl_ccb.dev.supercube2015.typ_a, 70
 hvl_ccb.dev.technix, 173
 hvl_ccb.dev.tiepie, 83
 hvl_ccb.dev.tiepie.base, 72
 hvl_ccb.dev.tiepie.channel, 74
 hvl_ccb.dev.tiepie.device, 77
 hvl_ccb.dev.tiepie.generator, 78
 hvl_ccb.dev.tiepie.i2c, 79
 hvl_ccb.dev.tiepie.oscilloscope, 80
 hvl_ccb.dev.tiepie.utils, 83
 hvl_ccb.dev.utils, 177
 hvl_ccb.dev.visa, 179
 hvl_ccb.experiment_manager, 185
 hvl_ccb.utils, 183
 hvl_ccb.utils.enum, 181
 hvl_ccb.utils.typing, 182
 hvl_ccb.utils.validation, 182
 MONITOR_I (*FuGProbusVRegisterGroups* attribute), 108
 MONITOR_V (*FuGProbusVRegisterGroups* attribute), 108
 most_recent_error (*FuGProbusVConfigRegisters* property), 107
 motion_distance_per_full_step (*NewportSMC100PPConfig* attribute), 143
 motor_config (*NewportSMC100PPConfig* property), 143
 move_to_absolute_position() (*NewportSMC100PP* method), 139
 move_to_relative_position() (*NewportSMC100PP* method), 139
 move_wait_sec (*NewportSMC100PPConfig* attribute), 143
 MOVING (*NewportSMC100PP.StateMessages* attribute), 136
 MOVING (*NewportStates* attribute), 147
 MS_NOMINAL_CURRENT (*PSI9000* attribute), 96
 MS_NOMINAL_VOLTAGE (*PSI9000* attribute), 96
 msb_first() (*TechnixStatusByte* method), 176
 MULTI_COMMANDS_MAX (*VisaCommunication* attribute), 30
 MULTI_COMMANDS_SEPARATOR (*VisaCommunication* attribute), 30

N

n_channels (*TiePieOscilloscope* property), 81
 n_max_try_get_device (*TiePieDeviceConfig* attribute), 73
 NameEnum (*class* in *hvl_ccb.utils.enum*), 181
 NAMES (*SerialCommunicationParity* attribute), 25
 names() (*RTO1024.TriggerModes* class method), 155
 namespace_index (*SupercubeConfiguration* attribute), 37, 60
 NATIVEONLY (*TiePieOscilloscopeAutoResolutionModes* attribute), 82
 NED_END_OF_TURN (*NewportSMC100PP.MotorErrors* attribute), 135
 NEG (*ILS2T.Ref16Jog* attribute), 163
 NEG_FAST (*ILS2T.Ref16Jog* attribute), 163
 NEGATIVE (*FuGPolarities* attribute), 105
 negative_software_limit (*NewportSMC100PPConfig* attribute), 143
 NewportConfigCommands (*class* in *hvl_ccb.dev.newport*), 134
 NewportControllerError, 135
 NewportMotorError, 135
 NewportMotorPowerSupplyWasCutError, 135
 NewportSerialCommunicationError, 147
 NewportSMC100PP (*class* in *hvl_ccb.dev.newport*), 135
 NewportSMC100PP.MotorErrors (*class* in *hvl_ccb.dev.newport*), 135
 NewportSMC100PP.StateMessages (*class* in *hvl_ccb.dev.newport*), 136
 NewportSMC100PPConfig (*class* in *hvl_ccb.dev.newport*), 141
 NewportSMC100PPConfig.EspStageConfig (*class* in *hvl_ccb.dev.newport*), 142
 NewportSMC100PPConfig.HomeSearch (*class* in *hvl_ccb.dev.newport*), 142
 NewportSMC100PPSerialCommunication (*class* in *hvl_ccb.dev.newport*), 143
 NewportSMC100PPSerialCommunication.ControllerErrors (*class* in *hvl_ccb.dev.newport*), 144
 NewportSMC100PPSerialCommunicationConfig (*class* in *hvl_ccb.dev.newport*), 146
 NewportStates (*class* in *hvl_ccb.dev.newport*), 147
 NewportUncertainPositionError, 147

NO (*FuGDigitalVal* attribute), 103
 NO_ERROR (*NewportSMC100PPSerialCommunication*.*Control* attribute), 144
 NO_REF (*NewportStates* attribute), 147
 NO_REF_ESP_STAGE_ERROR (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_CONFIG (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_DISABLED (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_HOMING (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_JOGGING (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_MOVING (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_READY (*portSMC100PP.StateMessages* attribute), 136
 NO_REF_FROM_RESET (*portSMC100PP.StateMessages* attribute), 136
 No_sensor (*PfeifferTPG.SensorStatus* attribute), 148
 NOISE (*TiePieGeneratorSignalType* attribute), 79
 NONE (*ILS2T.Ref16Jog* attribute), 163
 NONE (*LabJack.ThermocoupleType* attribute), 120
 None (*PfeifferTPG.SensorTypes* attribute), 149
 NONE (*SerialCommunicationParity* attribute), 25
 NoPower (*PowerSetup* attribute), 51
 NORMAL (*RTO1024.TriggerModes* attribute), 155
 noSen (*PfeifferTPG.SensorTypes* attribute), 149
 noSENSOR (*PfeifferTPG.SensorTypes* attribute), 149
 not_defined (*AlarmText* attribute), 40, 63
 nr_trials_activate (*LuminoxConfig* attribute), 170
 NullCommunicationProtocol (class in *hvl_ccb.comm.base*), 14
 number (*EarthingStick* property), 46
 Number (in module *hvl_ccb.utils.typing*), 182
 number_of_decimals (*HeinzingerConfig* attribute), 113
 number_of_sensors (*PfeifferTPG* property), 150

O

ODD (*SerialCommunicationParity* attribute), 25
 OFF (*FuGDigitalVal* attribute), 103
 OFF (*HeinzingerDI.OutputStatus* attribute), 113
 offset (*TiePieGeneratorConfig* property), 78
 OH (*NewportConfigCommands* attribute), 135
 Ok (*PfeifferTPG.SensorStatus* attribute), 149

on (*FuG property*), 101
 ON (*FuG* attribute), 103
 on (*FuGProbusVDIRegisters* property), 107
 ON (*HeinzingerDI.OutputStatus* attribute), 113
 ON (*ILS2T.State* attribute), 163
 ONE (*HeinzingerConfig.RecordingsEnum* attribute), 112
 ONE (*LabJack.AInRange* attribute), 118
 ONE (*SerialCommunicationStopbits* attribute), 25
 ONE_HUNDREDTH (*LabJack.AInRange* attribute), 118
 ONE_POINT_FIVE (*SerialCommunicationStopbits* attribute), 25
 ONE_TENTH (*LabJack.AInRange* attribute), 119
 ONLYUPWARDSOFTZERO (*FuGRampModes* attribute), 109
 OpcControl (class in *hvl_ccb.dev.supercube.constants*), 50
 OpcUaCommunication (class in *hvl_ccb.comm.opc*), 20
 OpcUaCommunicationConfig (class in *hvl_ccb.comm.opc*), 21
 OpcUaCommunicationIOError, 22
 OpcUaCommunicationTimeoutError, 22
 OpcUaSubHandler (class in *hvl_ccb.comm.opc*), 22
 open (*DoorStatus* attribute), 45, 63
 open (*EarthingStickOperation* attribute), 47
 open (*EarthingStickStatus* attribute), 47, 65
 open() (*CommunicationProtocol* method), 14
 open() (*LaudaProRp245eTcpCommunication* method), 129
 open() (*LJMCommunication* method), 15
 open() (*ModbusTcpCommunication* method), 18
 open() (*NullCommunicationProtocol* method), 14
 open() (*OpcUaCommunication* method), 20
 open() (*SerialCommunication* method), 23
 open() (*Tcp* method), 26
 open() (*TelnetCommunication* method), 28
 open() (*VisaCommunication* method), 30
 open_shutter() (*CryLasLaser* method), 91
 open_timeout (*VisaCommunicationConfig* attribute), 32
 OPENED (*CryLasLaser.AnswersShutter* attribute), 90
 OPENED (*CryLasLaserShutterStatus* attribute), 96
 operate() (*Supercube2015Base* method), 59
 operate() (*SupercubeBase* method), 35
 operate_earthing_stick() (*SupercubeBase* method), 35
 operating_status() (*EarthingStick* class method), 46
 operating_status_1 (*EarthingStick* attribute), 46
 operating_status_2 (*EarthingStick* attribute), 46
 operating_status_3 (*EarthingStick* attribute), 46
 operating_status_4 (*EarthingStick* attribute), 46
 operating_status_5 (*EarthingStick* attribute), 46
 operating_status_6 (*EarthingStick* attribute), 46
 operating_statuses() (*EarthingStick* class method), 46

OPERATION_MODE (*LaudaProRp245eCommand* attribute), 127
operation_mode (*LaudaProRp245eConfig* attribute), 129
optional_defaults() (*AsyncCommunicationProtocolConfig* class method), 13
optional_defaults() (*CryLasAttenuatorConfig* class method), 88
optional_defaults() (*CryLasAttenuatorSerialCommunicationConfig* class method), 89
optional_defaults() (*CryLasLaserConfig* class method), 93
optional_defaults() (*CryLasLaserSerialCommunicationConfig* class method), 95
optional_defaults() (*EmptyConfig* class method), 86, 184
optional_defaults() (*FuGConfig* class method), 103
optional_defaults() (*FuGSerialCommunicationConfig* class method), 111
optional_defaults() (*HeinzingerConfig* class method), 113
optional_defaults() (*HeinzingerSerialCommunicationConfig* class method), 117
optional_defaults() (*ILS2TConfig* class method), 166
optional_defaults() (*ILS2TModbusTcpCommunicationConfig* class method), 167
optional_defaults() (*LaudaProRp245eConfig* class method), 129
optional_defaults() (*LaudaProRp245eTcpCommunicationConfig* class method), 130
optional_defaults() (*LJMCommunicationConfig* class method), 17
optional_defaults() (*LuminoxConfig* class method), 170
optional_defaults() (*LuminoxSerialCommunicationConfig* class method), 172
optional_defaults() (*MBW973Config* class method), 132
optional_defaults() (*MBW973SerialCommunicationConfig* class method), 134
optional_defaults() (*ModbusTcpCommunicationConfig* class method), 19
optional_defaults() (*NewportSMC100PPConfig* class method), 143
optional_defaults() (*NewportSMC100PPSerialCommunicationConfig* class method), 147
optional_defaults() (*OpcUaCommunicationConfig* class method), 21
optional_defaults() (*PfeifferTPGConfig* class method), 151
optional_defaults() (*PfeifferTPGSerialCommunicationConfig* class method), 153
optional_defaults() (*PSI9000Config* class method), 99
optional_defaults() (*PSI9000VisaCommunicationConfig* class method), 100
optional_defaults() (*RTO1024Config* class method), 160
optional_defaults() (*RTO1024VisaCommunicationConfig* class method), 161
optional_defaults() (*SerialCommunicationConfig* class method), 25
optional_defaults() (*SupercubeAOpcUaConfiguration* class method), 54, 71
optional_defaults() (*SupercubeBOpcUaConfiguration* class method), 56
optional_defaults() (*SupercubeConfiguration* class method), 37, 60
optional_defaults() (*SupercubeOpcUaCommunicationConfig* class method), 38, 61
optional_defaults() (*TcpCommunicationConfig* class method), 27
optional_defaults() (*TechnixCommunicationConfig* class method), 174
optional_defaults() (*TechnixConfig* class method), 175
optional_defaults() (*TechnixSerialCommunicationConfig* class method), 176
optional_defaults() (*TechnixTelnetCommunicationConfig* class method), 177
optional_defaults() (*TelnetCommunicationConfig* class method), 29
optional_defaults() (*TiePieDeviceConfig* class method), 73
optional_defaults() (*VisaCommunicationConfig* class method), 32
optional_defaults() (*VisaDeviceConfig* class method), 180
OSCILLOSCOPE (*TiePieDeviceType* attribute), 73
OT (*NewportConfigCommands* attribute), 135
out (*FuGProbusVDORegisters* property), 108
out_1_1 (*GeneralSupport* attribute), 49, 66
out_1_2 (*GeneralSupport* attribute), 49, 66
out_2_1 (*GeneralSupport* attribute), 49, 66
out_2_2 (*GeneralSupport* attribute), 49, 66
out_3_1 (*GeneralSupport* attribute), 49, 66
out_3_2 (*GeneralSupport* attribute), 49, 66
out_4_1 (*GeneralSupport* attribute), 49, 66
out_4_2 (*GeneralSupport* attribute), 49, 66
out_5_1 (*GeneralSupport* attribute), 49, 66
out_5_2 (*GeneralSupport* attribute), 49, 66

out_6_1 (*GeneralSupport attribute*), 49, 66
out_6_2 (*GeneralSupport attribute*), 49, 66
OUTPUT (*FuGProbusIVCommands attribute*), 106
output() (*GeneralSupport class method*), 49, 66
output_off() (*FuGProbusIV method*), 105
output_off() (*HeinzingerDI method*), 114
output_on() (*HeinzingerDI method*), 114
OUTPUT_POWER_EXCEEDED (*New-portSMC100PP.MotorErrors attribute*), 135
output_status (*HeinzingerDI property*), 114
OUTPUTONCMD (*FuGProbusVRegisterGroups attribute*), 108
OUTPUTX0 (*FuGProbusVRegisterGroups attribute*), 108
OUTPUTX1 (*FuGProbusVRegisterGroups attribute*), 108
OUTPUTX2 (*FuGProbusVRegisterGroups attribute*), 108
OUTPUTXCMD (*FuGProbusVRegisterGroups attribute*), 109
outX0 (*FuG property*), 102
outX1 (*FuG property*), 102
outX2 (*FuG property*), 102
outXCMD (*FuG property*), 102
Overrange (*PfeifferTPG.SensorStatus attribute*), 149

P

PARAM_MISSING_OR_INVALID (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
parity (*CryLasAttenuatorSerialCommunicationConfig attribute*), 89
parity (*CryLasLaserSerialCommunicationConfig attribute*), 95
parity (*FuGSerialCommunicationConfig attribute*), 111
parity (*HeinzingerSerialCommunicationConfig attribute*), 117
parity (*LuminoxSerialCommunicationConfig attribute*), 172
parity (*MBW973SerialCommunicationConfig attribute*), 134
parity (*NewportSMC100PPSerialCommunicationConfig attribute*), 147
parity (*PfeifferTPGSerialCommunicationConfig attribute*), 153
Parity (*SerialCommunicationConfig attribute*), 24
parity (*SerialCommunicationConfig attribute*), 25
parse_read_measurement_value() (*LuminoxMeasurementType method*), 170
partial_pressure_o2 (*LuminoxMeasurementType attribute*), 170
Pascal (*PfeifferTPG.PressureUnits attribute*), 148
pause() (*LaudaProRp245e method*), 125
pause_ramp() (*LaudaProRp245e method*), 125
PBR (*PfeifferTPG.SensorTypes attribute*), 149
PEAK_CURRENT_LIMIT (*New-portSMC100PP.MotorErrors attribute*), 135
peak_output_current_limit (*New-portSMC100PPConfig attribute*), 143
percent_o2 (*LuminoxMeasurementType attribute*), 170
PfeifferTPG (*class in hvl_ccb.dev.pfeiffer_tpg*), 148
PfeifferTPG.PressureUnits (*class in hvl_ccb.dev.pfeiffer_tpg*), 148
PfeifferTPG.SensorStatus (*class in hvl_ccb.dev.pfeiffer_tpg*), 148
PfeifferTPG.SensorTypes (*class in hvl_ccb.dev.pfeiffer_tpg*), 149
PfeifferTPGConfig (*class in hvl_ccb.dev.pfeiffer_tpg*), 151
PfeifferTPGConfig.Model (*class in hvl_ccb.dev.pfeiffer_tpg*), 151
PfeifferTPGError, 152
PfeifferTPGSerialCommunication (*class in hvl_ccb.dev.pfeiffer_tpg*), 152
PfeifferTPGSerialCommunicationConfig (*class in hvl_ccb.dev.pfeiffer_tpg*), 152
PKR (*PfeifferTPG.SensorTypes attribute*), 149
POLARITY (*FuGProbusIVCommands attribute*), 106
Poller (*class in hvl_ccb.dev.utils*), 177
polling (*LuminoxOutputMode attribute*), 171
polling_delay_sec (*SupercubeConfiguration attribute*), 37
polling_interval (*MBW973Config attribute*), 133
polling_interval_sec (*SupercubeConfiguration attribute*), 37
polling_interval_sec (*TechnixConfig attribute*), 175
polling_period (*CryLasLaserConfig attribute*), 93
polling_timeout (*CryLasLaserConfig attribute*), 93
port (*ModbusTcpCommunicationConfig attribute*), 19
port (*OpcUaCommunicationConfig attribute*), 22
port (*SerialCommunicationConfig attribute*), 25
port (*SupercubeOpcUaCommunicationConfig attribute*), 61
port (*TcpCommunicationConfig attribute*), 27
port (*TechnixTelnetCommunicationConfig attribute*), 177
port (*TelnetCommunicationConfig attribute*), 29
port (*VisaCommunicationConfig attribute*), 32
port_range() (*GeneralSupport class method*), 49
POS (*ILS2T.Ref16Jog attribute*), 163
POS_END_OF_TURN (*NewportSMC100PP.MotorErrors attribute*), 135
POS_FAST (*ILS2T.Ref16Jog attribute*), 163
POSITION (*ILS2TRegAddr attribute*), 167
POSITION_OUT_OF_LIMIT (*New-portSMC100PPSerialCommunication.ControllerErrors attribute*), 144
POSITIVE (*FuGPolarities attribute*), 105

<code>positive_software_limit</code>	(<i>NewportSMC100PPConfig attribute</i>), 143	<code>query_multiple()</code>	(<i>NewportSMC100PPSerialCommunication method</i>), 145
<code>post_force_value()</code>	(<i>NewportSMC100PPConfig method</i>), 143	<code>query_polling()</code> (<i>Luminox method</i>), 169	
<code>post_stop_pause_sec</code> (<i>TechnixConfig attribute</i>), 175		<code>QUICKSTOP</code> (<i>ILS2T.State attribute</i>), 163	
<code>Power</code> (<i>class in hvl_ccb.dev.supercube.constants</i>), 50		<code>QuickStop</code> (<i>SafetyStatus attribute</i>), 52, 69	
<code>Power</code> (<i>class in hvl_ccb.dev.supercube2015.constants</i>), 67		<code>quickstop()</code> (<i>ILS2T method</i>), 164	
<code>power_limit</code> (<i>PSI9000Config attribute</i>), 99		<code>quit</code> (<i>Errors attribute</i>), 48, 65	
<code>PowerSetup</code> (<i>class in hvl_ccb.dev.supercube.constants</i>), 51		<code>quit_error()</code> (<i>Supercube2015Base method</i>), 59	
<code>PowerSetup</code> (<i>class in hvl_ccb.dev.supercube2015.constants</i>), 67		<code>quit_error()</code> (<i>SupercubeBase method</i>), 35	
<code>pre_sample_ratio</code> (<i>TiePieOscilloscopeConfig property</i>), 82			
<code>prepare_ultra_segmentation()</code> (<i>RTO1024 method</i>), 157			
<code>probe_offset</code> (<i>TiePieOscilloscopeChannelConfig property</i>), 75			
<code>PSI9000</code> (<i>class in hvl_ccb.dev.ea_psi9000</i>), 96			
<code>PSI9000Config</code> (<i>class in hvl_ccb.dev.ea_psi9000</i>), 98			
<code>PSI9000Error</code> , 99			
<code>PSI9000VisaCommunication</code> (<i>class in hvl_ccb.dev.ea_psi9000</i>), 99			
<code>PSI9000VisaCommunicationConfig</code> (<i>class in hvl_ccb.dev.ea_psi9000</i>), 99			
<code>PT100</code> (<i>LabJack.ThermocoupleType attribute</i>), 120			
<code>PT1000</code> (<i>LabJack.ThermocoupleType attribute</i>), 120			
<code>PT500</code> (<i>LabJack.ThermocoupleType attribute</i>), 120			
<code>PTP</code> (<i>ILS2T.Mode attribute</i>), 162			
<code>PublicPropertiesReprMixin</code> (<i>class in hvl_ccb.dev.tiepie.utils</i>), 83			
<code>PULSE</code> (<i>TiePieGeneratorSignalType attribute</i>), 79			
<code>pump_init</code> (<i>LaudaProRp245eConfig attribute</i>), 129			
<code>PUMP_LEVEL</code> (<i>LaudaProRp245eCommand attribute</i>), 127			
Q			
<code>QIL</code> (<i>NewportConfigCommands attribute</i>), 135			
<code>QUERY</code> (<i>FuGProbusIVCommands attribute</i>), 106			
<code>query()</code> (<i>CryLasLaserSerialCommunication method</i>), 94			
<code>query()</code> (<i>FuGSerialCommunication method</i>), 110			
<code>query()</code> (<i>NewportSMC100PPSerialCommunication method</i>), 145			
<code>query()</code> (<i>PfeifferTPGSerialCommunication method</i>), 152			
<code>query()</code> (<i>SyncCommunicationProtocol method</i>), 14			
<code>query()</code> (<i>TechnixCommunication method</i>), 174			
<code>query()</code> (<i>VisaCommunication method</i>), 30			
<code>query_all()</code> (<i>CryLasLaserSerialCommunication method</i>), 94			
<code>query_command()</code>	(<i>LaudaProRp245eTcpCommunication method</i>), 129		
		<code>read()</code> (<i>MBW973 method</i>), 131	
		<code>read()</code> (<i>OpcUaCommunication method</i>), 20	
		<code>read()</code> (<i>Supercube2015Base method</i>), 59	
		<code>read()</code> (<i>SupercubeBase method</i>), 35	
		<code>read()</code> (<i>Tcp method</i>), 26	
		<code>read_all()</code> (<i>AsyncCommunicationProtocol method</i>), 12	
		<code>read_bytes()</code> (<i>AsyncCommunicationProtocol method</i>), 12	
		<code>read_bytes()</code> (<i>SerialCommunication method</i>), 23	
		<code>read_bytes()</code> (<i>TelnetCommunication method</i>), 28	
		<code>read_float()</code> (<i>MBW973 method</i>), 131	
		<code>read_holding_registers()</code> (<i>ModbusTcpCommunication method</i>), 18	

```

read_input_registers() (ModbusTcpCommunication method), 18
read_int() (MBW973 method), 131
read_measurement() (RTO1024 method), 157
read_measurements() (MBW973 method), 131
read_name() (LJMCommunication method), 15
read_nonempty() (AsyncCommunicationProtocol method), 12
read_resistance() (LabJack method), 122
read_single_bytes() (SerialCommunication method), 23
read_streaming() (Luminox method), 169
read_termination (VisaCommunicationConfig attribute), 32
read_text() (AsyncCommunicationProtocol method), 12
read_text() (NewportSMC100PPSerialCommunication method), 145
read_text_nonempty() (AsyncCommunicationProtocol method), 12
READ_TEXT_SKIP_PREFIXES (CryLasLaserSerialCommunication attribute), 94
read_thermocouple() (LabJack method), 122
readback_data (FuGProbusVConfigRegisters property), 107
READBACKCHANNEL (FuGProbusIVCommands attribute), 106
READY (CryLasLaser.AnswersStatus attribute), 90
READY (ILS2T.State attribute), 163
READY (NewportStates attribute), 147
ready() (Supercube2015Base method), 59
ready() (SupercubeBase method), 35
READY_ACTIVE (CryLasLaser.LaserStatus attribute), 90
READY_FROM_DISABLE (New-portSMC100PP.StateMessages attribute), 136
READY_FROM_HOMING (New-portSMC100PP.StateMessages attribute), 136
READY_FROM_JOGGING (New-portSMC100PP.StateMessages attribute), 136
READY_FROM_MOVING (New-portSMC100PP.StateMessages attribute), 136
READY_INACTIVE (CryLasLaser.LaserStatus attribute), 90
record_length (TiePieOscilloscopeConfig property), 82
RedOperate (SafetyStatus attribute), 52, 69
RedReady (SafetyStatus attribute), 52, 69
reg_3 (FuGProbusVDIRegisters property), 108
RegAddr (ILS2T attribute), 163
RegDatatype (ILS2T attribute), 163
register_pulse_time (TechnixConfig attribute), 175
RELATIVE (TiePieOscilloscopeTriggerLevelMode attribute), 76
RELATIVE_POSITION_MOTOR (ILS2T.ActionsPtp attribute), 162
RELATIVE_POSITION_TARGET (ILS2T.ActionsPtp attribute), 162
remote (Technix property), 173
remove_device() (DeviceSequenceMixin method), 85
require_block_measurement_support (TiePieDeviceConfig attribute), 73
required_keys() (AsyncCommunicationProtocolConfig class method), 13
required_keys() (CryLasAttenuatorConfig class method), 88
required_keys() (CryLasAttenuatorSerialCommunicationConfig class method), 89
required_keys() (CryLasLaserConfig class method), 93
required_keys() (CryLasLaserSerialCommunicationConfig class method), 95
required_keys() (EmptyConfig class method), 86, 184
required_keys() (FuGConfig class method), 103
required_keys() (FuGSerialCommunicationConfig class method), 111
required_keys() (HeinzingerConfig class method), 113
required_keys() (HeinzingerSerialCommunicationConfig class method), 117
required_keys() (ILS2TConfig class method), 166
required_keys() (ILS2TModbusTcpCommunicationConfig class method), 167
required_keys() (LaudaProRp245eConfig class method), 129
required_keys() (LaudaProRp245eTcpCommunicationConfig class method), 130
required_keys() (LJMCommunicationConfig class method), 17
required_keys() (LuminoxConfig class method), 170
required_keys() (LuminoxSerialCommunicationConfig class method), 172
required_keys() (MBW973Config class method), 133
required_keys() (MBW973SerialCommunicationConfig class method), 134
required_keys() (ModbusTcpCommunicationConfig class method), 19
required_keys() (NewportSMC100PPConfig class method), 143
required_keys() (New-portSMC100PPSerialCommunicationConfig class method), 147
required_keys() (OpcUaCommunicationConfig class method), 22

```

required_keys() (*PfeifferTPGConfig class method*), 151
required_keys() (*PfeifferTPGSerialCommunicationConfig class method*), 153
required_keys() (*PSI9000Config class method*), 99
required_keys() (*PSI9000VisaCommunicationConfig class method*), 100
required_keys() (*RTO1024Config class method*), 161
required_keys() (*RTO1024VisaCommunicationConfig class method*), 161
required_keys() (*SerialCommunicationConfig class method*), 25
required_keys() (*SupercubeAOpcUaConfiguration class method*), 54, 72
required_keys() (*SupercubeBOpcUaConfiguration class method*), 56
required_keys() (*SupercubeConfiguration class method*), 37, 61
required_keys() (*SupercubeOpcUaCommunicationConfig class method*), 38, 61
required_keys() (*TcpCommunicationConfig class method*), 27
required_keys() (*TechnixCommunicationConfig class method*), 174
required_keys() (*TechnixConfig class method*), 175
required_keys() (*TechnixSerialCommunicationConfig class method*), 176
required_keys() (*TechnixTelnetCommunicationConfig class method*), 177
required_keys() (*TelnetCommunicationConfig class method*), 29
required_keys() (*TiePieDeviceConfig class method*), 73
required_keys() (*VisaCommunicationConfig class method*), 32
required_keys() (*VisaDeviceConfig class method*), 180
reset (*BreakdownDetection attribute*), 45, 63
RESET (*FuGProbusIVCommands attribute*), 106
reset() (*FuGProbusIV method*), 106
reset() (*NewportSMC100PP method*), 140
reset() (*VisaDevice method*), 179
reset_error() (*ILS2T method*), 164
reset_interface() (*HeinzingerDI method*), 114
reset_ramp() (*LaudaProRp245e method*), 125
resolution (*TiePieOscilloscopeConfig property*), 82
response_sleep_time (*CryLasAttenuatorConfig attribute*), 88
RISING (*TiePieOscilloscopeTriggerKind attribute*), 76
RISING_OR_FALLING (*TiePieOscilloscopeTriggerKind attribute*), 76
RMS_CURRENT_LIMIT (*NewportSMC100PP.MotorErrors attribute*), 135
rpm_max_init (*ILS2TConfig attribute*), 166
rs485_address (*NewportSMC100PPConfig attribute*), 143
RT01024 (*class in hvl_ccb.dev.rs_rto1024*), 154
RT01024.TriggerModes (*class in hvl_ccb.dev.rs_rto1024*), 154
RT01024Config (*class in hvl_ccb.dev.rs_rto1024*), 160
RT01024Error, 161
RT01024VisaCommunication (*class in hvl_ccb.dev.rs_rto1024*), 161
RT01024VisaCommunicationConfig (*class in hvl_ccb.dev.rs_rto1024*), 161
run() (*ExperimentManager method*), 186
run() (*LaudaProRp245e method*), 125
run_continuous_acquisition() (*RTO1024 method*), 157
run_single_acquisition() (*RTO1024 method*), 157
RUNNING (*ExperimentStatus attribute*), 186

S

S (*LabJack.ThermocoupleType attribute*), 120
SA (*NewportConfigCommands attribute*), 135
safe_ground_enabled (*TiePieOscilloscopeChannelConfig property*), 75
Safety (*class in hvl_ccb.dev.supercube.constants*), 51
Safety (*class in hvl_ccb.dev.supercube2015.constants*), 68
SafetyStatus (*class in hvl_ccb.dev.supercube.constants*), 52
SafetyStatus (*class in hvl_ccb.dev.supercube2015.constants*), 68
sample_frequency (*TiePieOscilloscopeConfig property*), 82
save_configuration() (*RTO1024 method*), 157
save_waveform_history() (*RTO1024 method*), 157
SCALE (*ILS2TRegAddr attribute*), 167
ScalingFactorValueError, 168
screw_scaling (*NewportSMC100PPConfig attribute*), 143
send_command() (*NewportSMC100PPSerialCommunication method*), 145
send_command() (*PfeifferTPGSerialCommunication method*), 152
send_pulses() (*LabJack method*), 122
send_stop() (*NewportSMC100PPSerialCommunication method*), 146
Sensor_error (*PfeifferTPG.SensorStatus attribute*), 149
Sensor_off (*PfeifferTPG.SensorStatus attribute*), 149
sensor_status (*LuminoxMeasurementType attribute*), 171
SERIAL (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 128
serial_number (*LuminoxMeasurementType attribute*), 171

<code>serial_number (TiePieDeviceConfig attribute)</code> , 73			<code>set_motor_configuration()</code> (<i>NewportSMC100PP method</i>), 140
<code>SerialCommunication (class in hvl_ccb.comm.serial)</code> , 23			<code>set_negative_software_limit()</code> (<i>NewportSMC100PP method</i>), 140
<code>SerialCommunicationBytesize</code> (class) <i>in</i> <code>hvl_ccb.comm.serial</code> , 24			<code>set_number_of_recordings()</code> (<i>HeinzingerDI method</i>), 114
<code>SerialCommunicationConfig</code> (class) <i>in</i> <code>hvl_ccb.comm.serial</code> , 24			<code>set_output()</code> (<i>PSI9000 method</i>), 97
<code>SerialCommunicationIOError</code> , 25			<code>set_positive_software_limit()</code> (<i>NewportSMC100PP method</i>), 140
<code>SerialCommunicationParity</code> (class) <i>in</i> <code>hvl_ccb.comm.serial</code> , 25			<code>set_pulse_energy()</code> (<i>CryLasLaser method</i>), 91
<code>SerialCommunicationStopbits</code> (class) <i>in</i> <code>hvl_ccb.comm.serial</code> , 25			<code>set_pump_level()</code> (<i>LaudaProRp245e method</i>), 125
<code>set_acceleration()</code> (<i>NewportSMC100PP method</i>), 140			<code>set_ramp_iterations()</code> (<i>LaudaProRp245e method</i>), 125
<code>set_acquire_length()</code> (<i>RTO1024 method</i>), 158			<code>set_ramp_program()</code> (<i>LaudaProRp245e method</i>), 126
<code>set_ain_differential()</code> (<i>LabJack method</i>), 122			<code>set_ramp_segment()</code> (<i>LaudaProRp245e method</i>), 126
<code>set_ain_range()</code> (<i>LabJack method</i>), 122			<code>set_ramp_type()</code> (<i>ILS2T method</i>), 165
<code>set_ain_resistance()</code> (<i>LabJack method</i>), 123			<code>set_reference_point()</code> (<i>RTO1024 method</i>), 159
<code>set_ain_resolution()</code> (<i>LabJack method</i>), 123			<code>set_register()</code> (<i>FuGProbusV method</i>), 106
<code>set_ain_thermocouple()</code> (<i>LabJack method</i>), 123			<code>set_remote_control()</code> (<i>Supercube2015Base method</i>), 59
<code>set_analog_output()</code> (<i>LabJack method</i>), 123			<code>set_remote_control()</code> (<i>SupercubeBase method</i>), 36
<code>set_attenuation()</code> (<i>CryLasAttenuator method</i>), 87			<code>set_repetition_rate()</code> (<i>CryLasLaser method</i>), 92
<code>set_cee16_socket()</code> (<i>Supercube2015Base method</i>), 59			<code>set_repetitions()</code> (<i>RTO1024 method</i>), 159
<code>set_cee16_socket()</code> (<i>SupercubeBase method</i>), 35			<code>set_slope()</code> (<i>Supercube2015WithFU method</i>), 71
<code>set_channel_offset()</code> (<i>RTO1024 method</i>), 158			<code>set_slope()</code> (<i>SupercubeWithFU method</i>), 55
<code>set_channel_position()</code> (<i>RTO1024 method</i>), 158			<code>set_status_board()</code> (<i>SupercubeBase method</i>), 36
<code>set_channel_range()</code> (<i>RTO1024 method</i>), 158			<code>set_support_output()</code> (<i>Supercube2015Base method</i>), 59
<code>set_channel_scale()</code> (<i>RTO1024 method</i>), 158			<code>set_support_output()</code> (<i>SupercubeBase method</i>), 36
<code>set_channel_state()</code> (<i>RTO1024 method</i>), 159			<code>set_support_output_impulse()</code> (<i>Supercube2015Base method</i>), 59
<code>set_clock()</code> (<i>LabJack method</i>), 123			<code>set_support_output_impulse()</code> (<i>SupercubeBase method</i>), 36
<code>set_control_mode()</code> (<i>LaudaProRp245e method</i>), 125			<code>set_system_lock()</code> (<i>PSI9000 method</i>), 97
<code>set_current()</code> (<i>HeinzingerDI method</i>), 114			<code>set_t13_socket()</code> (<i>Supercube2015Base method</i>), 60
<code>set_current()</code> (<i>HeinzingerPNC method</i>), 115			<code>set_t13_socket()</code> (<i>SupercubeBase method</i>), 36
<code>set_digital_output()</code> (<i>LabJack method</i>), 124			<code>set_target_voltage()</code> (<i>Supercube2015WithFU method</i>), 71
<code>set_display_unit()</code> (<i>PfeifferTPG method</i>), 150			<code>set_target_voltage()</code> (<i>SupercubeWithFU method</i>), 55
<code>set_earthing_manual()</code> (<i>Supercube2015Base method</i>), 59			<code>set_temp_set_point()</code> (<i>LaudaProRp245e method</i>), 126
<code>set_external_temp()</code> (<i>LaudaProRp245e method</i>), 125			<code>set_transmission()</code> (<i>CryLasAttenuator method</i>), 87
<code>set_full_scale_mbar()</code> (<i>PfeifferTPG method</i>), 150			<code>set_trigger_level()</code> (<i>RTO1024 method</i>), 159
<code>set_full_scale_unitless()</code> (<i>PfeifferTPG method</i>), 150			<code>set_trigger_mode()</code> (<i>RTO1024 method</i>), 160
<code>set_init_attenuation()</code> (<i>CryLasAttenuator method</i>), 87			<code>set_trigger_source()</code> (<i>RTO1024 method</i>), 160
<code>set_init_shutter_status()</code> (<i>CryLasLaser method</i>), 91			<code>set_upper_limits()</code> (<i>PSI9000 method</i>), 98
<code>set_jog_speed()</code> (<i>ILS2T method</i>), 164			<code>set_voltage()</code> (<i>HeinzingerDI method</i>), 114
<code>set_lower_limits()</code> (<i>PSI9000 method</i>), 97			<code>set_voltage()</code> (<i>HeinzingerPNC method</i>), 115
<code>set_max_acceleration()</code> (<i>ILS2T method</i>), 165			<code>set_voltage_current()</code> (<i>PSI9000 method</i>), 98
<code>set_max_deceleration()</code> (<i>ILS2T method</i>), 165			<code>SETCURRENT</code> (<i>FuGProbusVRegisterGroups attribute</i>), 109
<code>set_max_rpm()</code> (<i>ILS2T method</i>), 165			<code>setup</code> (<i>Power attribute</i>), 51, 67
<code>set_measuring_options()</code> (<i>MBW973 method</i>), 132			<code>setvalue</code> (<i>FuGProbusVSetRegisters property</i>), 109
<code>set_message_board()</code> (<i>SupercubeBase method</i>), 36			

SETVOLTAGE (*FuGProbusVRegisterGroups attribute*), 109
 SEVENBITS (*SerialCommunicationBytesize attribute*), 24
 SHORT_CIRCUIT (*NewportSMC100PP.MotorErrors attribute*), 136
 SHUTDOWN_CURRENT_LIMIT (*PSI9000 attribute*), 96
 SHUTDOWN_VOLTAGE_LIMIT (*PSI9000 attribute*), 96
 ShutterStatus (*CryLasLaser attribute*), 90
 ShutterStatus (*CryLasLaserConfig attribute*), 93
 signal_type (*TiePieGeneratorConfig property*), 78
 SINE (*TiePieGeneratorSignalType attribute*), 79
 SingleCommDevice (*class in hvl_ccb.dev.base*), 86
 SIXBITS (*SerialCommunicationBytesize attribute*), 24
 SIXTEEN (*HeinzingerConfig.RecordingsEnum attribute*), 112
 SIXTEEN_BIT (*TiePieOscilloscopeResolution attribute*), 82
 SL (*NewportConfigCommands attribute*), 135
 SN (*FuGReadbackChannels attribute*), 110
 SOFTWARE_INTERNAL_SIXTY (*CryLasLaser.RepetitionRates attribute*), 90
 SOFTWARE_INTERNAL_TEN (*CryLasLaser.RepetitionRates attribute*), 90
 SOFTWARE_INTERNAL_TWENTY (*CryLasLaser.RepetitionRates attribute*), 90
 software_revision (*LuminoxMeasurementType attribute*), 171
 SPACE (*SerialCommunicationParity attribute*), 25
 SPECIALRAMPUPWARDS (*FuGRampModes attribute*), 109
 spoll() (*VisaCommunication method*), 31
 spoll_handler() (*VisaDevice method*), 179
 SQUARE (*TiePieGeneratorSignalType attribute*), 79
 SR (*NewportConfigCommands attribute*), 135
 srq_mask (*FuGProbusVConfigRegisters property*), 107
 srq_status (*FuGProbusVConfigRegisters property*), 107
 stage_configuration (*NewportSMC100PPConfig attribute*), 143
 START (*LaudaProRp245eCommand attribute*), 127
 start() (*CryLasAttenuator method*), 87
 start() (*CryLasLaser method*), 92
 start() (*Device method*), 84
 start() (*DeviceSequenceMixin method*), 85
 start() (*ExperimentManager method*), 186
 start() (*FuG method*), 102
 start() (*FuGProbusIV method*), 106
 start() (*HeinzingerDI method*), 114
 start() (*HeinzingerPNC method*), 115
 start() (*ILS2T method*), 165
 start() (*LabJack method*), 124
 start() (*LaudaProRp245e method*), 126
 start() (*Luminox method*), 169
 start() (*MBW973 method*), 132
 start() (*NewportSMC100PP method*), 141
 start() (*PfeifferTPG method*), 150
 start() (*PSI9000 method*), 98
 start() (*RTO1024 method*), 160
 start() (*SingleCommDevice method*), 86
 start() (*Supercube2015Base method*), 60
 start() (*SupercubeBase method*), 37
 start() (*Technix method*), 173
 start() (*TiePieGeneratorMixin method*), 79
 start() (*TiePieI2CHostMixin method*), 80
 start() (*TiePieOscilloscope method*), 81
 start() (*VisaDevice method*), 180
 start_control() (*MBW973 method*), 132
 start_measurement() (*TiePieOscilloscope method*), 81
 start_polling() (*Poller method*), 178
 start_ramp() (*LaudaProRp245e method*), 126
 STARTING (*ExperimentStatus attribute*), 186
 States (*NewportSMC100PP attribute*), 136
 status (*ExperimentManager property*), 186
 status (*FuGProbusVConfigRegisters property*), 107
 status (*FuGProbusVDORegisters property*), 108
 status (*Safety attribute*), 52
 status() (*EarthingStick class method*), 46
 status_1 (*Door attribute*), 45
 status_1 (*EarthingRod attribute*), 45
 status_1 (*EarthingStick attribute*), 47
 status_1_closed (*EarthingStick attribute*), 64
 status_1_connected (*EarthingStick attribute*), 64
 status_1_open (*EarthingStick attribute*), 64
 status_2 (*Door attribute*), 45
 status_2 (*EarthingRod attribute*), 45
 status_2 (*EarthingStick attribute*), 47
 status_2_closed (*EarthingStick attribute*), 64
 status_2_connected (*EarthingStick attribute*), 64
 status_2_open (*EarthingStick attribute*), 64
 status_3 (*Door attribute*), 45
 status_3 (*EarthingRod attribute*), 45
 status_3 (*EarthingStick attribute*), 47
 status_3_closed (*EarthingStick attribute*), 64
 status_3_connected (*EarthingStick attribute*), 64
 status_3_open (*EarthingStick attribute*), 64
 status_4 (*EarthingStick attribute*), 47
 status_4_closed (*EarthingStick attribute*), 64
 status_4_connected (*EarthingStick attribute*), 64
 status_4_open (*EarthingStick attribute*), 64
 status_5 (*EarthingStick attribute*), 47
 status_5_closed (*EarthingStick attribute*), 64
 status_5_connected (*EarthingStick attribute*), 64
 status_5_open (*EarthingStick attribute*), 64
 status_6 (*EarthingStick attribute*), 47
 status_6_closed (*EarthingStick attribute*), 64
 status_6_connected (*EarthingStick attribute*), 64
 status_6_open (*EarthingStick attribute*), 64
 status_closed() (*EarthingStick class method*), 64

status_connected() (*EarthingStick class method*), 64
status_error (*Safety attribute*), 68
status_green (*Safety attribute*), 68
status_open() (*EarthingStick class method*), 65
status_ready_for_red (*Safety attribute*), 68
status_red (*Safety attribute*), 68
STATUSBYTE (*FuGReadbackChannels attribute*), 110
statuses() (*EarthingStick class method*), 47
stop (*Errors attribute*), 48, 65
STOP (*LaudaProRp245eCommand attribute*), 127
stop() (*CryLasLaser method*), 92
stop() (*Device method*), 84
stop() (*DeviceSequenceMixin method*), 85
stop() (*ExperimentManager method*), 186
stop() (*FuGProbusIV method*), 106
stop() (*HeinzingerDI method*), 115
stop() (*ILS2T method*), 165
stop() (*LabJack method*), 124
stop() (*LaudaProRp245e method*), 126
stop() (*Luminox method*), 169
stop() (*MBW973 method*), 132
stop() (*NewportSMC100PP method*), 141
stop() (*PfeifferTPG method*), 151
stop() (*PSI9000 method*), 98
stop() (*RTO1024 method*), 160
stop() (*SingleCommDevice method*), 86
stop() (*Supercube2015Base method*), 60
stop() (*SupercubeBase method*), 37
stop() (*Technix method*), 174
stop() (*TiePieGeneratorMixin method*), 79
stop() (*TiePieI2CHostMixin method*), 80
stop() (*TiePieOscilloscope method*), 81
stop() (*VisaDevice method*), 180
stop_acquisition() (*RTO1024 method*), 160
stop_motion() (*NewportSMC100PP method*), 141
stop_number (*Errors attribute*), 65
stop_polling() (*Poller method*), 178
stop_ramp() (*LaudaProRp245e method*), 126
stopbits (*CryLasAttenuatorSerialCommunicationConfig attribute*), 89
stopbits (*CryLasLaserSerialCommunicationConfig attribute*), 95
stopbits (*FuGSerialCommunicationConfig attribute*), 111
stopbits (*HeinzingerSerialCommunicationConfig attribute*), 117
stopbits (*LuminoxSerialCommunicationConfig attribute*), 172
stopbits (*MBW973SerialCommunicationConfig attribute*), 134
stopbits (*NewportSMC100PPSerialCommunicationConfig attribute*), 147
stopbits (*PfeifferTPGSerialCommunicationConfig attribute*), 153
Stopbits (*SerialCommunicationConfig attribute*), 24
stopbits (*SerialCommunicationConfig attribute*), 25
streaming (*LuminoxOutputMode attribute*), 171
StrEnumBase (*class in hvl_ccb.utils.enum*), 181
sub_handler (*OpcUaCommunicationConfig attribute*), 22
sub_handler (*SupercubeOpcUaCommunicationConfig attribute*), 38, 61
suitable_range() (*TiePieOscilloscopeRange static method*), 76
Supercube2015Base (*class in hvl_ccb.dev.supercube2015.base*), 57
Supercube2015WithFU (*class in hvl_ccb.dev.supercube2015.typ_a*), 70
SupercubeA0pcUaCommunication (*class in hvl_ccb.dev.supercube.typ_a*), 53
SupercubeA0pcUaCommunication (*class in hvl_ccb.dev.supercube2015.typ_a*), 71
SupercubeA0pcUaConfiguration (*class in hvl_ccb.dev.supercube.typ_a*), 53
SupercubeA0pcUaConfiguration (*class in hvl_ccb.dev.supercube2015.typ_a*), 71
SupercubeB (*class in hvl_ccb.dev.supercube.typ_b*), 55
SupercubeBase (*class in hvl_ccb.dev.supercube.base*), 33
SupercubeB0pcUaCommunication (*class in hvl_ccb.dev.supercube.typ_b*), 56
SupercubeB0pcUaConfiguration (*class in hvl_ccb.dev.supercube.typ_b*), 56
SupercubeConfiguration (*class in hvl_ccb.dev.supercube.base*), 37
SupercubeConfiguration (*class in hvl_ccb.dev.supercube2015.base*), 60
SupercubeEarthlingStickOperationError, 38
SupercubeOpcEndpoint (*class in hvl_ccb.dev.supercube.constants*), 52
SupercubeOpcEndpoint (*class in hvl_ccb.dev.supercube2015.constants*), 69
SupercubeOpcUaCommunication (*class in hvl_ccb.dev.supercube.base*), 38
SupercubeOpcUaCommunication (*class in hvl_ccb.dev.supercube2015.base*), 61
SupercubeOpcUaCommunicationConfig (*class in hvl_ccb.dev.supercube.base*), 38
SupercubeOpcUaCommunicationConfig (*class in hvl_ccb.dev.supercube2015.base*), 61
SupercubeSubscriptionHandler (*class in hvl_ccb.dev.supercube.base*), 38
SupercubeSubscriptionHandler (*class in hvl_ccb.dev.supercube2015.base*), 62
SupercubeWithFU (*class in hvl_ccb.dev.supercube.typ_a*), 54
switch_to_operate (*Safety attribute*), 52
switch_to_ready (*Safety attribute*), 52

<code>switchto_green</code> (<i>Safety attribute</i>), 68					
<code>switchto_operate</code> (<i>Safety attribute</i>), 68					
<code>switchto_ready</code> (<i>Safety attribute</i>), 68					
<code>SyncCommunicationProtocol</code> (class in <i>hvl_ccb.comm.base</i>), 14					
<code>SyncCommunicationProtocolConfig</code> (class in <i>hvl_ccb.comm.base</i>), 14					
T					
<code>T</code> (<i>LabJack.ThermocoupleType attribute</i>), 120					
<code>t13_1</code> (<i>GeneralSockets attribute</i>), 48, 65					
<code>t13_2</code> (<i>GeneralSockets attribute</i>), 48, 65					
<code>t13_3</code> (<i>GeneralSockets attribute</i>), 48, 65					
<code>T13_SOCKET_PORTS</code> (in module <i>hvl_ccb.dev.supercube.constants</i>), 52					
<code>T13_SOCKET_PORTS</code> (in module <i>hvl_ccb.dev.supercube2015.constants</i>), 69					
<code>T1MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T200MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T20MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T256US</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T4</code> (<i>LabJack.DeviceType attribute</i>), 120					
<code>T4 (LJMCommunicationConfig.DeviceType attribute)</code> , 16					
<code>T40MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T4MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T7</code> (<i>LabJack.DeviceType attribute</i>), 120					
<code>T7 (LJMCommunicationConfig.DeviceType attribute)</code> , 16					
<code>T7_PRO</code> (<i>LabJack.DeviceType attribute</i>), 120					
<code>T7_PRO</code> (<i>LJMCommunicationConfig.DeviceType attribute</i>), 16					
<code>T800MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>T80MS</code> (<i>FuGMonitorModes attribute</i>), 105					
<code>target_pulse_energy</code> (<i>CryLasLaser property</i>), 92					
<code>Tcp</code> (class in <i>hvl_ccb.comm.tcp</i>), 26					
<code>TCP</code> (<i>LJMCommunicationConfig.ConnectionType attribute</i>), 16					
<code>TcpCommunicationConfig</code> (class in <i>hvl_ccb.comm.tcp</i>), 26					
<code>TCP/IP_INSTR</code> (<i>VisaCommunicationConfig.InterfaceType attribute</i>), 31					
<code>TCPIP_SOCKET</code> (<i>VisaCommunicationConfig.InterfaceType attribute</i>), 31					
<code>TEC1</code> (<i>CryLasLaser.AnswersStatus attribute</i>), 90					
<code>TEC2</code> (<i>CryLasLaser.AnswersStatus attribute</i>), 90					
<code>Technix</code> (class in <i>hvl_ccb.dev.technix</i>), 173					
<code>TechnixCommunication</code> (class in <i>hvl_ccb.dev.technix</i>), 174					
<code>TechnixCommunicationConfig</code> (class in <i>hvl_ccb.dev.technix</i>), 174					
<code>TechnixConfig</code> (class in <i>hvl_ccb.dev.technix</i>), 174					
<code>TechnixError</code> , 175					
<code>TechnixSerialCommunication</code> (class in <i>hvl_ccb.dev.technix</i>), 175					
			<code>TechnixSerialCommunicationConfig</code> (class in <i>hvl_ccb.dev.technix</i>), 176		
			<code>TechnixStatusByte</code> (class in <i>hvl_ccb.dev.technix</i>), 176		
			<code>TechnixTelnetCommunication</code> (class in <i>hvl_ccb.dev.technix</i>), 176		
			<code>TechnixTelnetCommunicationConfig</code> (class in <i>hvl_ccb.dev.technix</i>), 177		
			<code>TelnetCommunication</code> (class in <i>hvl_ccb.comm.telnet</i>), 28		
			<code>TelnetCommunicationConfig</code> (class in <i>hvl_ccb.comm.telnet</i>), 29		
			<code>TelnetError</code> , 29		
			<code>TEMP</code> (<i>ILS2TRegAddr attribute</i>), 167		
			<code>TEMP_SET_POINT</code> (<i>LaudaProRp245eCommand attribute</i>), 127		
			<code>temp_set_point_init</code> (<i>LaudaProRp245eConfig attribute</i>), 129		
			<code>temperature_sensor</code> (<i>LuminoxMeasurementType attribute</i>), 171		
			<code>TEN</code> (<i>LabJack.AInRange attribute</i>), 119		
			<code>TEN</code> (<i>LabJack.CalMicroAmpere attribute</i>), 119		
			<code>TEN_MHZ</code> (<i>LabJack.ClockFrequency attribute</i>), 119		
			<code>terminator</code> (<i>AsyncCommunicationProtocolConfig attribute</i>), 14		
			<code>terminator</code> (<i>CryLasAttenuatorSerialCommunicationConfig attribute</i>), 89		
			<code>terminator</code> (<i>CryLasLaserSerialCommunicationConfig attribute</i>), 96		
			<code>TERMINATOR</code> (<i>FuGProbusIVCommands attribute</i>), 106		
			<code>terminator</code> (<i>FuGProbusVConfigRegisters property</i>), 107		
			<code>terminator</code> (<i>FuGSerialCommunicationConfig attribute</i>), 111		
			<code>terminator</code> (<i>HeinzingerSerialCommunicationConfig attribute</i>), 117		
			<code>terminator</code> (<i>LaudaProRp245eTcpCommunicationConfig attribute</i>), 130		
			<code>terminator</code> (<i>LuminoxSerialCommunicationConfig attribute</i>), 172		
			<code>terminator</code> (<i>MBW973SerialCommunicationConfig attribute</i>), 134		
			<code>terminator</code> (<i>NewportSMC100PPSerialCommunicationConfig attribute</i>), 147		
			<code>terminator</code> (<i>PfeifferTPGSerialCommunicationConfig attribute</i>), 153		
			<code>terminator</code> (<i>TechnixCommunicationConfig attribute</i>), 174		
			<code>terminator_str()</code> (<i>SerialCommunicationConfig method</i>), 25		
			<code>THIRTY_TWO_BIT</code> (<i>LabJack.BitLimit attribute</i>), 119		
			<code>TiePieDeviceConfig</code> (class in <i>hvl_ccb.dev.tiepie.base</i>), 72		
			<code>TiePieDeviceType</code> (class in <i>hvl_ccb.dev.tiepie.base</i>), 73		

TiePieError	73		
TiePieGeneratorConfig	(class <i>hvl_ccb.dev.tiepie.generator</i>), 78	in	timeout (PfeifferTPGSerialCommunicationConfig attribute), 153
TiePieGeneratorConfigLimits	(class <i>hvl_ccb.dev.tiepie.generator</i>), 78	in	timeout (SerialCommunicationConfig attribute), 25
TiePieGeneratorMixin	(class <i>hvl_ccb.dev.tiepie.generator</i>), 78	in	timeout (TelnetCommunicationConfig attribute), 29
TiePieGeneratorSignalType	(class <i>hvl_ccb.dev.tiepie.generator</i>), 79	in	timeout (VisaCommunicationConfig attribute), 32
TiePieHS5	(class in <i>hvl_ccb.dev.tiepie.device</i>), 77		Torr (PfeifferTPG.PressureUnits attribute), 148
TiePieHS6	(class in <i>hvl_ccb.dev.tiepie.device</i>), 77		TPG25xA (PfeifferTPGConfig.Model attribute), 151
TiePieI2CHostConfig	(class in <i>hvl_ccb.dev.tiepie.i2c</i>), 79		TPGx6x (PfeifferTPGConfig.Model attribute), 151
TiePieI2CHostConfigLimits	(class <i>hvl_ccb.dev.tiepie.i2c</i>), 79	in	TPR (PfeifferTPG.SensorTypes attribute), 149
TiePieI2CHostMixin	(class in <i>hvl_ccb.dev.tiepie.i2c</i>), 79		transmission (CryLasAttenuator property), 87
TiePieOscilloscope	(class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 80	in	TRIANGLE (TiePieGeneratorSignalType attribute), 79
TiePieOscilloscopeAutoResolutionModes	(class in <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 81		trigger_enabled (TiePieOscilloscopeChannelConfig property), 75
TiePieOscilloscopeChannelConfig	(class in <i>hvl_ccb.dev.tiepie.channel</i>), 74	in	trigger_hysteresis (TiePieOscilloscopeChannel- Config property), 75
TiePieOscilloscopeChannelConfigLimits	(class in <i>hvl_ccb.dev.tiepie.channel</i>), 75		trigger_kind (TiePieOscilloscopeChannelConfig prop- erty), 75
TiePieOscilloscopeChannelCoupling	(class in <i>hvl_ccb.dev.tiepie.channel</i>), 75	in	trigger_level (TiePieOscilloscopeChannelConfig property), 75
TiePieOscilloscopeConfig	(class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 82	in	trigger_level_mode (TiePieOscilloscopeChannel- Config property), 75
TiePieOscilloscopeConfigLimits	(class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 82	in	trigger_timeout (TiePieOscilloscopeConfig prop- erty), 82
TiePieOscilloscopeRange	(class <i>hvl_ccb.dev.tiepie.channel</i>), 75	in	triggered (BreakdownDetection attribute), 45, 63
TiePieOscilloscopeResolution	(class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 82	in	TWELVE_BIT (TiePieOscilloscopeResolution attribute), 82
TiePieOscilloscopeTriggerKind	(class <i>hvl_ccb.dev.tiepie.channel</i>), 76	in	TWELVE_HUNDRED_FIFTY_KHZ (Lab- Jack.ClockFrequency attribute), 119
TiePieOscilloscopeTriggerLevelMode	(class in <i>hvl_ccb.dev.tiepie.channel</i>), 76		TWENTY_FIVE_HUNDRED_KHZ (LabJack.ClockFrequency attribute), 119
TiePieWS5	(class in <i>hvl_ccb.dev.tiepie.device</i>), 77		TWENTY_MHZ (LabJack.ClockFrequency attribute), 119
timeout (CryLasAttenuatorSerialCommunicationConfig attribute)	, 89		TWENTY_VOLT (TiePieOscilloscopeRange attribute), 76
timeout (CryLasLaserSerialCommunicationConfig at- tribute)	, 96		TWO (HeinzingerConfig.RecordingsEnum attribute), 112
timeout (FuGSerialCommunicationConfig attribute)	, 111		TWO (SerialCommunicationStopbits attribute), 26
timeout (HeinzingerSerialCommunicationConfig attribute)	, 117		TWO_HUNDRED (LabJack.CalMicroAmpere attribute), 119
timeout (LuminoxSerialCommunicationConfig at- tribute)	, 172		TWO_HUNDRED_MILLI_VOLT (TiePieOscilloscopeRange attribute), 76
timeout (MBW973SerialCommunicationConfig at- tribute)	, 134		TWO_VOLT (TiePieOscilloscopeRange attribute), 76
timeout (NewportSMC100PPSerialCommunicationConfig attribute)	, 147		U
			Underrange (PfeifferTPG.SensorStatus attribute), 149
			unit (ILS2TModbusTcpCommunicationConfig attribute), 167
			unit (ModbusTcpCommunicationConfig attribute), 19
			unit (PfeifferTPG property), 151
			unit_current (HeinzingerPNC property), 115
			unit_voltage (HeinzingerPNC property), 115
			UNKNOWN (HeinzingerDI.OutputStatus attribute), 113
			UNKNOWN (HeinzingerPNC.UnitCurrent attribute), 115
			UNKNOWN (HeinzingerPNC.UnitVoltage attribute), 115
			UNKNOWN (TiePieGeneratorSignalType attribute), 79
			UNKNOWN (TiePieOscilloscopeAutoResolutionModes at- tribute), 82

UNKNOWN (*TiePieOscilloscopeTriggerLevelMode attribute*), 76

UNREADY_INACTIVE (*CryLasLaser.LaserStatus attribute*), 90

update_laser_status() (*CryLasLaser method*), 92

update_period (*OpcUaCommunicationConfig attribute*), 22

update_repetition_rate() (*CryLasLaser method*), 92

update_shutter_status() (*CryLasLaser method*), 92

update_target_pulse_energy() (*CryLasLaser method*), 92

UpdateEspStageInfo (*NewportSMC100PPConfig.EspStageConfig attribute*), 142

UPPER_TEMP (*LaudaProRp245eCommand attribute*), 127

upper_temp (*LaudaProRp245eConfig attribute*), 129

USB (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 128

USB (*LJMCommunicationConfig.ConnectionType attribute*), 16

user_position_offset (*NewportSMC100PPConfig attribute*), 143

user_steps() (*ILS2T method*), 165

V

V (*HeinzingerPNC.UnitVoltage attribute*), 115

VA (*NewportConfigCommands attribute*), 135

validate_bool() (*in module hvl_ccb.utils.validation*), 182

validate_number() (*in module hvl_ccb.utils.validation*), 182

validate_pump_level() (*LaudaProRp245e method*), 126

value (*FuGProbusVMonitorRegisters property*), 108

value (*LabJack.AInRange property*), 119

value_raw (*FuGProbusVMonitorRegisters property*), 108

ValueEnum (*class in hvl_ccb.utils.enum*), 181

VB (*NewportConfigCommands attribute*), 135

velocity (*NewportSMC100PPConfig attribute*), 143

visa_backend (*VisaCommunicationConfig attribute*), 32

VisaCommunication (*class in hvl_ccb.comm.visa*), 30

VisaCommunicationConfig (*class in hvl_ccb.comm.visa*), 31

VisaCommunicationConfig.InterfaceType (*class in hvl_ccb.comm.visa*), 31

VisaCommunicationError, 32

VisaDevice (*class in hvl_ccb.dev.visa*), 179

VisaDeviceConfig (*class in hvl_ccb.dev.visa*), 180

VOLT (*ILS2TRegAddr attribute*), 167

Volt (*PfeifferTPG.PressureUnits attribute*), 148

voltage (*FuG property*), 102

VOLTAGE (*FuGProbusIVCommands attribute*), 106

VOLTAGE (*FuGReadbackChannels attribute*), 110

voltage (*Technix property*), 174

voltage_lower_limit (*PSI9000Config attribute*), 99

voltage_max (*Power attribute*), 51, 67

voltage_monitor (*FuG property*), 102

voltage_primary (*Power attribute*), 51, 67

voltage_regulation (*Technix property*), 174

voltage_slope (*Power attribute*), 51, 67

voltage_target (*Power attribute*), 51, 67

voltage_upper_limit (*PSI9000Config attribute*), 99

W

WAIT_AFTER_WRITE (*VisaCommunication attribute*), 30

wait_for_polling_result() (*Poller method*), 178

wait_operation_complete() (*VisaDevice method*), 180

wait_sec_initialisation (*PSI9000Config attribute*), 99

wait_sec_max_disable (*ILS2TConfig attribute*), 166

wait_sec_post_absolute_position (*ILS2TConfig attribute*), 166

wait_sec_post_activate (*LuminoxConfig attribute*), 170

wait_sec_post_CANNOT_DISABLE (*ILS2TConfig attribute*), 166

wait_sec_post_enable (*ILS2TConfig attribute*), 166

wait_sec_post_relative_step (*ILS2TConfig attribute*), 166

wait_sec_pre_read_or_write (*LaudaProRp245eTcpCommunicationConfig attribute*), 130

wait_sec_read_text_nonempty (*AsyncCommunicationProtocolConfig attribute*), 14

wait_sec_read_text_nonempty (*FuGSerialCommunicationConfig attribute*), 111

wait_sec_read_text_nonempty (*HeinzingerSerialCommunicationConfig attribute*), 117

wait_sec_retry_get_device (*TiePieDeviceConfig attribute*), 73

wait_sec_settings_effect (*PSI9000Config attribute*), 99

wait_sec_stop_commands (*FuGConfig attribute*), 103

wait_sec_stop_commands (*HeinzingerConfig attribute*), 113

wait_sec_system_lock (*PSI9000Config attribute*), 99

wait_sec_trials_activate (*LuminoxConfig attribute*), 170

wait_timeout_retry_sec (*OpcUaCommunicationConfig attribute*), 22

wait_until_motor_initialized() (*NewportSMC100PP method*), 141

wait_until_ready() (*CryLasLaser method*), 93

warning (*Errors attribute*), 48

WIFI (*LJMCommunicationConfig.ConnectionType attribute*), 16
wrap_libtiepie_exception() (*in module hvl_ccb.dev.tiepie.base*), 74
write() (*AsyncCommunicationProtocol method*), 12
write() (*MBW973 method*), 132
write() (*OpcUaCommunication method*), 21
write() (*Supercube2015Base method*), 60
write() (*SupercubeBase method*), 37
write() (*Tcp method*), 26
write() (*VisaCommunication method*), 31
write_absolute_position() (*ILS2T method*), 165
write_bytes() (*AsyncCommunicationProtocol method*), 13
write_bytes() (*SerialCommunication method*), 23
write_bytes() (*TelnetCommunication method*), 28
write_command() (*LaudaProRp245eTcpCommunication method*), 130
write_name() (*LJMCommunication method*), 16
write_names() (*LJMCommunication method*), 16
write_registers() (*ModbusTcpCommunication method*), 18
write_relative_step() (*ILS2T method*), 165
write_termination (*VisaCommunicationConfig attribute*), 32
write_text() (*AsyncCommunicationProtocol method*), 13
WRONG_ESP_STAGE (*NewportSMC100PP.MotorErrors attribute*), 136

X

x_stat (*FuGProbusVDIRegisters property*), 108
XOUTPUTS (*FuGProbusIVCommands attribute*), 106

Y

YES (*FuGDigitalVal attribute*), 103

Z

ZX (*NewportConfigCommands attribute*), 135