
HVL Common Code Base Documentation

Release 0.9.0

Mikolaj Rybiński, David Graber, Henrik Menne, Alise Chachereau,

Jan 07, 2022

CONTENTS:

1	HVL Common Code Base	1
1.1	Features	1
1.2	Documentation	5
1.3	Credits	5
2	Installation	7
2.1	Stable release	7
2.2	From sources	7
2.3	Additional system libraries	8
3	Usage	9
4	API Documentation	11
4.1	hvl_ccb	11
5	Contributing	195
5.1	Types of Contributions	195
5.2	Get Started!	196
5.3	Merge Request Guidelines	197
5.4	Tips	197
5.5	Deploying	198
6	Credits	199
6.1	Maintainers	199
6.2	Authors	199
6.3	Contributors	199
7	History	201
7.1	0.9.0 (2022-01-07)	201
7.2	0.8.5 (2021-11-05)	201
7.3	0.8.4 (2021-10-22)	202
7.4	0.8.3 (2021-09-27)	202
7.5	0.8.2 (2021-08-27)	202
7.6	0.8.1 (2021-08-13)	202
7.7	0.8.0 (2021-07-02)	202
7.8	0.7.1 (2021-06-04)	203
7.9	0.7.0 (2021-05-25)	203
7.10	0.6.1 (2021-05-08)	203
7.11	0.6.0 (2021-04-23)	204
7.12	0.5.0 (2020-11-11)	204
7.13	0.4.0 (2020-07-16)	205

7.14	0.3.5 (2020-02-18)	205
7.15	0.3.4 (2019-12-20)	206
7.16	0.3.3 (2019-05-08)	206
7.17	0.3.2 (2019-05-08)	206
7.18	0.3.1 (2019-05-02)	206
7.19	0.3 (2019-05-02)	206
7.20	0.2.1 (2019-04-01)	207
7.21	0.2.0 (2019-03-31)	207
7.22	0.1.0 (2019-02-06)	207
8	Indices and tables	209
	Python Module Index	211
	Index	213

**CHAPTER
ONE**

HVL COMMON CODE BASE

Python common code base (CCB) to control devices, which are used in high-voltage research. All implemented devices are used and tested in the High Voltage Laboratory ([HVL](#)) of the Federal Institute of Technology Zurich (ETH Zurich).

- Free software: GNU General Public License v3
- Copyright (c) 2019-2022 ETH Zurich, SIS ID and HVL D-ITET

1.1 Features

For managing multi-device experiments instantiate the `ExperimentManager` utility class.

1.1.1 Devices

The device wrappers in `hvl_ccb` provide a standardised API with configuration dataclasses, various settings and options, as well as start/stop methods. Currently wrappers are available to control the following devices:

Function/Type	Devices
Data acquisition	LabJack (T4, T7, T7-PRO; requires LJM Library) Pico Technology PT-104 Platinum Resistance Data Logger (requires PicoSDK/libusbpt104)
Digital Delay Generator	Highland T560
Digital IO	LabJack (T4, T7, T7-PRO; requires LJM Library)
Experiment control	HVL Supercube with and without Frequency Converter
Gas Analyser	MBW 973-SF6 gas dew point mirror analyzer Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges SST Luminox oxygen sensor
I2C host	TiePie (HS5, WS5; requires LibTiePie SDK)
Laser	CryLaS pulsed laser CryLaS laser attenuator
Oscilloscope	Rhode & Schwarz RTO 1024 TiePie (HS5, HS6, WS5; requires LibTiePie SDK)
Power supply	Elektro-Automatik PSI9000 FuG Elektronik Heinzinger PNC Technix capacitor charger
Stepper motor drive	Newport SMC100PP Schneider Electric ILS2T
Temperature control	Lauda PRO RP 245 E circulation thermostat
Waveform generator	TiePie (HS5, WS5; requires LibTiePie SDK)
2	Chapter 1. HVL Common Code Base

Each device uses at least one standardised communication protocol wrapper.

1.1.2 Communication protocols

In hvl_ccb by “communication protocol” we mean different levels of communication standards, from the low level actual communication protocols like serial communication to application level interfaces like VISA TCP standard. There are also devices in hvl_ccb that use a dummy communication protocol; this is because these devices are build on proprietary manufacturer libraries that communicate with the corresponding devices, as in the case of TiePie or LabJack devices.

The communication protocol wrappers in hvl_ccb provide a standardised API with configuration dataclasses, as well as open/close and read/write/query methods. Currently, wrappers for the following communication protocols are available:

Communication protocol	Devices using
Modbus TCP	Schneider Electric ILS2T stepper motor drive
OPC UA	HVL Supercube with and without Frequency Converter
Serial	CryLaS pulsed laser and laser attenuator FuG Elektronik power supply (e.g. capacitor charger HCK) using the Probus V protocol Heinzinger PNC power supply using Heinzinger Digital Interface I/II SST Luminox oxygen sensor MBW 973-SF6 gas dew point mirror analyzer Newport SMC100PP single axis driver for 2-phase stepper motors Pfeiffer Vacuum TPG (25x, 26x and 36x) controller for compact pressure gauges Technix capacitor charger
TCP	Lauda PRO RP 245 E circulation thermostat
Telnet	Technix capacitor charger
VISA TCP	Elektro-Automatik PSI9000 DC power supply Rhode & Schwarz RTO 1024 oscilloscope
<i>proprietary</i>	LabJack (T4, T7, T7-PRO) devices, which communicate via LJM Library Pico Technology PT-104 Platinum Resistance Data Logger, which communicate via PicoSDK/libusbpt104 TiePie (HS5, HS6, WS5) oscilloscopes, generators and I2C hosts, which communicate via LibTiePie SDK

1.1.3 Sensor and Unit Conversion Utility

The Sensor and Unit Conversion Utility is a submodule that allows on the one hand a unified implementation of hardware-sensors and on the other hand provides a unified way to convert units. The utility can be used with single numbers (`int`, `float`) as well as array-like structures containing single numbers (`np.array()`, `list`, `dict`, `tuple`).

Currently the following sensors are implemented:

- LEM LT 4000S
- LMT 70A

The following unit conversion classes are implemented:

- Temperature (Kelvin, Celsius, Fahrenheit)
- Pressure (Pascal, Bar, Atmosphere, Psi, Torr, Millimeter Mercury)

1.2 Documentation

Note: if you're planning to contribute to the `hvl_ccb` project read the **Contributing** section in the HVL CCB documentation.

Do either:

- read [HVL CCB documentation at RTD](#),

or

- build and read HVL CCB documentation locally; install first [Graphviz](#) (make sure to have the `dot` command in the executable search path) and the Python build requirements for documentation:

```
$ pip install docs/requirements.txt
```

and then either on Windows in Git BASH run:

```
$ ./make.sh docs
```

or from any other shell with GNU Make installed run:

```
$ make docs
```

The target index HTML ("`docs/_build/html/index.html`") should open automatically in your Web browser.

1.3 Credits

This package was created with [Cookiecutter](#) and the [audreyr/cookiecutter-pypackage](#) project template.

INSTALLATION

2.1 Stable release

To install HVL Common Code Base, run this command in your terminal:

```
$ pip install hvl_ccb
```

To install HVL Common Code Base with optional Python libraries that require manual installations of additional system libraries, you need to specify on installation extra requirements corresponding to these controllers. For instance, to install Python requirements for LabJack and TiePie devices, run:

```
$ pip install "hvl_ccb[tiepie,labjack]"
```

See below for the info about additional system libraries and the corresponding extra requirements.

To install all extra requirements run:

```
$ pip install "hvl_ccb[all]"
```

This is the preferred method to install HVL Common Code Base, as it will always install the most recent stable release.

If you don't have `pip` installed, this [Python installation guide](#) can guide you through the process.

2.2 From sources

The sources for HVL Common Code Base can be downloaded from the [GitLab](#) repo.

You can either clone the repository:

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

Or download the [tarball](#):

```
$ curl -OL https://gitlab.com/ethz_hvl/hvl_ccb/-/archive/master/hvl_ccb.tar.gz
```

Once you have a copy of the source, you can install it with:

```
$ pip install .
```

2.3 Additional system libraries

If you have installed *hvl_ccb* with any of the extra features corresponding to device controllers, you must additionally install respective system library; these are:

Extra feature	Additional system library
labjack	LJM Library
picotech	PicoSDK (Windows) / libusbpt104 (Ubuntu/Debian)
tiepie	LibTiePie SDK

For more details on installation of the libraries see docstrings of the corresponding `hvl_ccb` modules.

**CHAPTER
THREE**

USAGE

To use HVL Common Code Base in a project:

```
import hvl_ccb
```


API DOCUMENTATION


4.1 hvl_ccb

4.1.1 Subpackages

[hvl_ccb.comm](#)

Submodules

[hvl_ccb.comm.base](#)

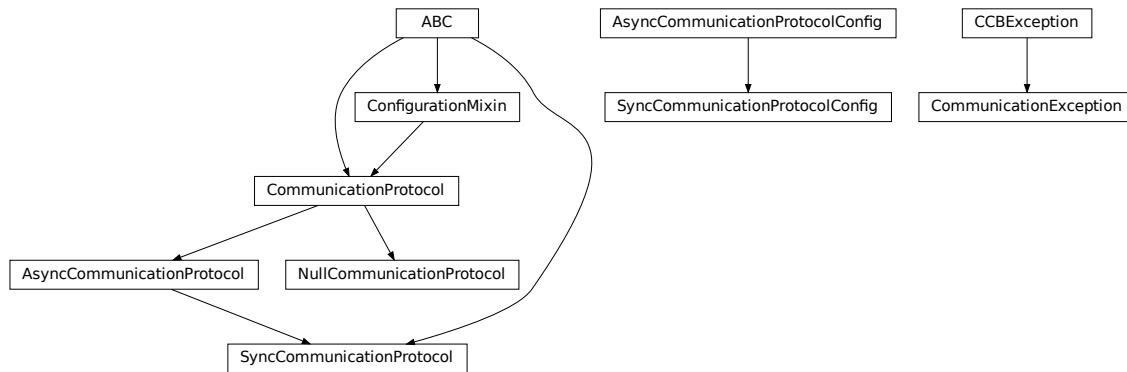

Module with base classes for communication protocols.

class `AsyncCommunicationProtocol(config)`

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Abstract base class for asynchronous communication protocols

static config_cls() → Type[`hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

read() → str

Read a single line of text as `str` from the communication.

Returns text as `str` including the terminator, which can also be empty “”

read_all(*n_attempts_max*: *Optional[int]* = *None*, *attempt_interval_sec*: *Optional[Union[int, float]]* = *None*) → *Optional[str]*

Read all lines of text from the connection till nothing is left to read.

Parameters

- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt_interval_sec** – time between the reading attempts

Returns A multi-line *str* including the terminator internally

abstract **read_bytes**() → *bytes*

Read a single line as *bytes* from the communication.

This method uses *self.access_lock* to ensure thread-safety.

Returns a single line as *bytes* containing the terminator, which can also be empty b””

read_nonempty(*n_attempts_max*: *Optional[int]* = *None*, *attempt_interval_sec*: *Optional[Union[int, float]]* = *None*) → *Optional[str]*

Try to read a non-empty single line of text as *str* from the communication. If the host does not reply or reply with white space only, it will return *None*.

Returns a non-empty text as a *str* or *None* in case of an empty string

Parameters

- **n_attempts_max** – Amount of attempts how often a non-empty text is tried to be read
- **attempt_interval_sec** – time between the reading attempts

read_text() → *str*

Read one line of text from the serial port. The input buffer may hold additional data afterwards, since only one line is read.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Returns String read from the serial port; “ if there was nothing to read.

Raises *SerialCommunicationIOError* – when communication port is not opened

read_text_nonempty(*n_attempts_max*: *Optional[int]* = *None*, *attempt_interval_sec*: *Optional[Union[int, float]]* = *None*) → *Optional[str]*

Reads from the serial port, until a non-empty line is found, or the number of attempts is exceeded.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Attention: in contrast to *read_text*, the returned answer will be stripped of a whitespace newline terminator at the end, if such terminator is set in the initial configuration (default).

Parameters

- **n_attempts_max** – maximum number of read attempts
- **attempt_interval_sec** – time between the reading attempts

Returns String read from the serial port; “ if number of attempts is exceeded or serial port is not opened.

write(*text*: *str*)

Write text as *str* to the communication.

Parameters **text** – test as a *str* to be written

abstract write_bytes(*data: bytes*) → int
 Write data as *bytes* to the communication.

This method uses *self.access_lock* to ensure thread-safety.

Parameters **data** – data as *bytes*-string to be written

Returns number of bytes written

write_text(*text: str*)
 Write text to the serial port. The text is encoded and terminated by the configured terminator.

NOTE: backward-compatibility proxy for *read* method; to be removed in v1.0

Parameters **text** – Text to send to the port.

Raises *SerialCommunicationIOError* – when communication port is not opened

class AsyncCommunicationProtocolConfig(*terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10*)

Bases: object

Base configuration data class for asynchronous communication protocols

clean_values()

default_n_attempts_read_text_nonempty: int = 10
 default number of attempts to read a non-empty text

encoding: str = 'utf-8'
 Standard encoding of the connection. Typically this is **utf-8**, but can also be **latin-1** or something from here: <https://docs.python.org/3/library/codecs.html#standard-encodings>

encoding_error_handling: str = 'strict'
 Encoding error handling scheme as defined here: <https://docs.python.org/3/library/codecs.html#error-handlers> By default strict error handling that raises *UnicodeError*.

force_value(*fieldname, value*)
 Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
 Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

terminator: bytes = b'\r\n'
The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations.
This defines the end of a single line.

wait_sec_read_text_nonempty: Union[int, float] = 0.5
time to wait between attempts of reading a non-empty text

exception CommunicationException

Bases: *hvl_ccb.exception.CCException*

class CommunicationProtocol(config)

Bases: *hvl_ccb.configuration.ConfigurationMixin, abc.ABC*

Communication protocol abstract base class.

Specifies the methods to implement for communication protocol, as well as implements some default settings and checks.

access_lock

Access lock to use with context manager when accessing the communication protocol (thread safety)

abstract close()

Close the communication protocol

abstract open()

Open communication protocol

class NullCommunicationProtocol(config)

Bases: *hvl_ccb.comm.base.CommunicationProtocol*

Communication protocol that does nothing.

close() → None

Void close function.

static config_cls() → Type[*hvl_ccb.configuration.EmptyConfig*]

Empty configuration

Returns EmptyConfig

open() → None

Void open function.

class SyncCommunicationProtocol(config)

Bases: *hvl_ccb.comm.base.AsyncCommunicationProtocol, abc.ABC*

Abstract base class for synchronous communication protocols with *query()*

static config_cls() → Type[*hvl_ccb.comm.base.SyncCommunicationProtocolConfig*]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(command: str) → Optional[str]


Send a command to the interface and handle the status message. Eventually raises an exception.

Parameters **command** – Command to send

Returns Answer from the interface, which can be None instead of an empty reply

```
class SyncCommunicationProtocolConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                      encoding_error_handling: str = 'strict',
                                      wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                      default_n_attempts_read_text_nonempty: int = 10)
Bases: hvl_ccb.comm.base.AsyncCommunicationProtocolConfig
```

hvl_ccb.comm.labjack_ljm

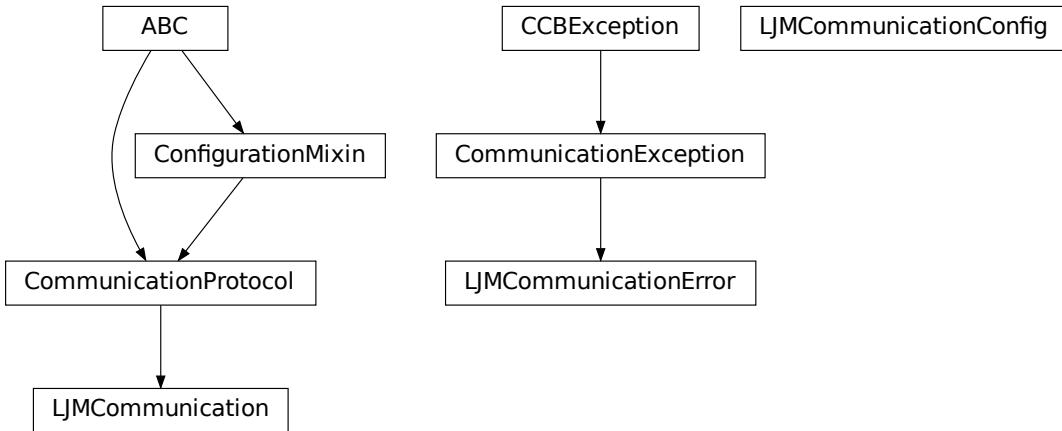

Communication protocol for LabJack using the LJM Library. Originally developed and tested for LabJack T7-PRO.

Makes use of the LabJack LJM Library Python wrapper. This wrapper needs an installation of the LJM Library for Windows, Mac OS X or Linux. Go to: <https://labjack.com/support/software/installers/ljm> and <https://labjack.com/support/software/examples/ljm/python>

```
class LJMCommunication(configuration)
Bases: hvl_ccb.comm.base.CommunicationProtocol
```

Communication protocol implementing the LabJack LJM Library Python wrapper.

close() → None
Close the communication port.

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

property is_open: bool
Flag indicating if the communication port is open.

Returns `True` if the port is open, otherwise `False`

open() → None
Open the communication port.

read_name(*names: str, return_num_type: Type[numbers.Real] = <class 'float'>) → Union[numbers.Real, Sequence[numbers.Real]]
Read one or more input numeric values by name.

Parameters

- **names** – one or more names to read out from the LabJack
- **return_num_type** – optional numeric type specification for return values; by default *float*.

Returns answer of the LabJack, either single number or multiple numbers in a sequence, respectively, when one or multiple names to read were given

Raises **TypeError** – if read value of type not compatible with *return_num_type*

write_name(*name*: str, *value*: numbers.Real) → None

Write one value to a named output.

Parameters

- **name** – String or with name of LabJack IO
- **value** – is the value to write to the named IO port

write_names(*name_value_dict*: Dict[str, numbers.Real]) → None

Write more than one value at once to named outputs.

Parameters **name_value_dict** – is a dictionary with string names of LabJack IO as keys and corresponding numeric values

```
class LJMCommunicationConfig(device_type: Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY',
                               connection_type: Union[str,
                               hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType] =
                               'ANY', identifier: str = 'ANY')
```

Bases: object

Configuration dataclass for [LJMCommunication](#).

```
class ConnectionType(value=<no_arg>, names=None, module=None, type=None, start=1,
                      boundary=None)
```

Bases: [hvl_ccb.utils.enum.AutoNumberNameEnum](#)

LabJack connection type.

ANY = 1

ETHERNET = 4

TCP = 3

USB = 2

WIFI = 5

```
class DeviceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: [hvl_ccb.utils.enum.AutoNumberNameEnum](#)

LabJack device types.

Can be also looked up by ambiguous Product ID (*p_id*) or by instance name: `python LabJackDeviceType(4)` is `LabJackDeviceType('T4')`

ANY = 1

T4 = 2

T7 = 3

T7_PRO = 4

classmethod `get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
List[hvl_ccb._dev.labjack.DeviceType]]`

Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.

Parameters `p_id` – Product ID of a LabJack device

Returns Instance or list of instances of `LabJackDeviceType`

Raises `ValueError` – when Product ID is unknown

clean_values() → None

Performs value checks on `device_type` and `connection_type`.

connection_type: `Union[str,`

`hvl_ccb.comm.labjack_ljm.LJMCommunicationConfig.ConnectionType] = 'ANY'`

Can be either string or of enum `ConnectionType`.

device_type: `Union[str, hvl_ccb._dev.labjack.DeviceType] = 'ANY'`

Can be either string 'ANY', 'T7_PRO', 'T7', 'T4', or of enum `DeviceType`.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

identifier: `str = 'ANY'`

The identifier specifies information for the connection to be used. This can be an IP address, serial number, or device name. See the LabJack docs (<https://labjack.com/support/software/api/ljm/function-reference/ljmopens/identifier-parameter>) for more information.

is_configdataclass = True

classmethod `keys()` → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod `optional_defaults()` → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod `required_keys()` → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns a list of strings containing all required keys.

exception `LJMCommunicationError`

Bases: `hvl_ccb.comm.base.CommunicationException`

Errors coming from LJMCommunication.

hvl_ccb.comm.modbus_tcp

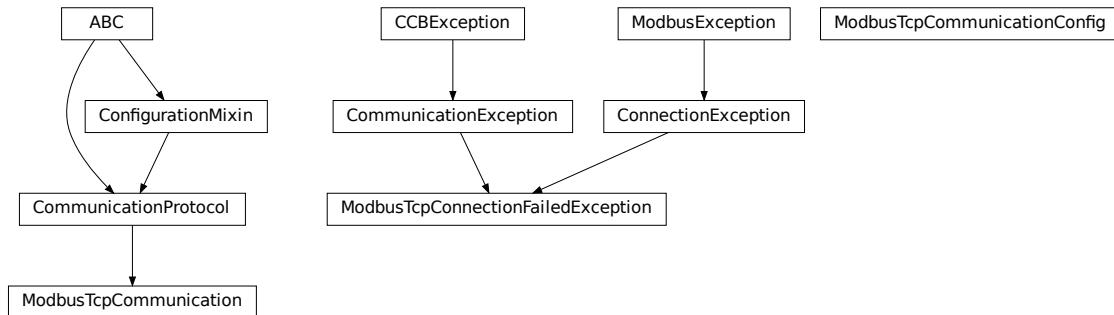

Communication protocol for modbus TCP ports. Makes use of the [pymodbus](#) library.

class **ModbusTcpCommunication**(*configuration*)
Bases: [hvl_ccb.comm.base.CommunicationProtocol](#)

Implements the Communication Protocol for modbus TCP.

close()

Close the Modbus TCP connection.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

open() → None

Open the Modbus TCP connection.

Raises [**ModbusTcpConnectionFailedException**](#) – if the connection fails.

read_holding_registers(*address*: int, *count*: int) → List[int]

Read specified number of register starting with given address and return the values from each register.

Parameters

- **address** – address of the first register
- **count** – count of registers to read

Returns list of int values

read_input_registers(*address*: int, *count*: int) → List[int]

Read specified number of register starting with given address and return the values from each register in a list.

Parameters

- **address** – address of the first register
- **count** – count of registers to read

Returns list of int values

write_registers(*address*: int, *values*: Union[List[int], int])

Write values from the specified address forward.

Parameters

- **address** – address of the first register
- **values** – list with all values

```
class ModbusTcpCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], unit: int, port: int = 502)
```

Bases: object

Configuration dataclass for [ModbusTcpCommunication](#).

clean_values()

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]
```

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 502

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

unit: int


```
exception ModbusTcpConnectionFailedException(string="")
```

Bases: pymodbus.exceptions.ConnectionException,

[hvl_ccb.comm.base.CommunicationException](#)

Exception raised when the connection failed.

`hvl_ccb.comm.opc`

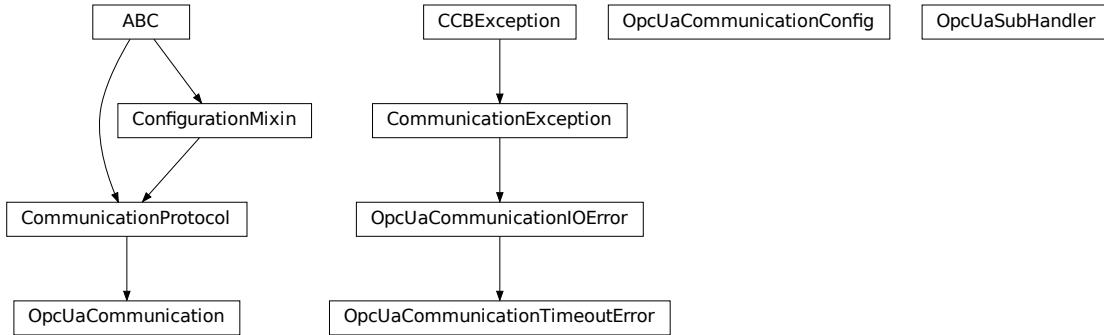

Communication protocol implementing an OPC UA connection. This protocol is used to interface with the “Super-cube” PLC from Siemens.

class `OpcUaCommunication(config)`

Bases: `hvl_ccb.comm.base.CommunicationProtocol`

Communication protocol implementing an OPC UA connection. Makes use of the package python-opcua.

`close() → None`

Close the connection to the OPC UA server.

`static config_cls()`

Return the default configdataclass class.

Returns a reference to the default configdataclass class

`init_monitored_nodes(node_id: Union[object, Iterable], ns_index: int) → None`

Initialize monitored nodes.

Parameters

- **`node_id`** – one or more strings of node IDs; node IDs are always casted via `str()` method here, hence do not have to be strictly string objects.
- **`ns_index`** – the namespace index the nodes belong to.

Raises `OpcUaCommunicationIOError` – when protocol was not opened or can't communicate with a OPC UA server

`property is_open: bool`

Flag indicating if the communication port is open.

Returns `True` if the port is open, otherwise `False`

`open() → None`

Open the communication to the OPC UA server.

Raises `OpcUaCommunicationIOError` – when communication port cannot be opened.

`read(node_id, ns_index)`

Read a value from a node with id and namespace index.

Parameters

- **`node_id`** – the ID of the node to read the value from

- **ns_index** – the namespace index of the node

Returns the value of the node object.

Raises *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

write(*node_id*, *ns_index*, *value*) → None

Write a value to a node with name *name*.

Parameters

- **node_id** – the id of the node to write the value to.
- **ns_index** – the namespace index of the node.
- **value** – the value to write.

Raises *OpcUaCommunicationIOError* – when protocol was not opened or can't communicate with a OPC UA server

```
class OpcUaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],  
                               endpoint_name: str, port: int = 4840, sub_handler:  
                               hvl_ccb.comm.opc.OpcUaSubHandler =  
                               <hvl_ccb.comm.opc.OpcUaSubHandler object>, update_period: int = 500,  
                               wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int =  
                               5)
```

Bases: *object*

Configuration dataclass for OPC UA Communciation.

clean_values()

endpoint_name: str

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]

Hostname or IP-Address of the OPC UA server.

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

max_timeout_retry_nr: int = 5

Maximal number of call re-tries on underlying OPC UA client timeout error

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

```

port: int = 4840
    Port of the OPC UA server to connect to.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

        Returns a list of strings containing all required keys.

sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.comm.opc.OpcUaSubHandler object>
    object to use for handling subscriptions.

update_period: int = 500
    Update period for generating datachange events in OPC UA [milli seconds]

wait_timeout_retry_sec: Union[int, float] = 1
    Wait time between re-trying calls on underlying OPC UA client timeout error

exception OpcUaCommunicationIOError
    Bases: OSError, hvl_ccb.comm.base.CommunicationException
    OPC-UA communication I/O error.

exception OpcUaCommunicationTimeoutError
    Bases: hvl_ccb.comm.opc.OpcUaCommunicationIOError
    OPC-UA communication timeout error.


class OpcUaSubHandler
    Bases: object
    Base class for subscription handling of OPC events and data change events. Override methods from this class to add own handling capabilities.

    To receive events from server for a subscription data_change and event methods are called directly from receiving thread. Do not do expensive, slow or network operation there. Create another thread if you need to do such a thing.

        datachange_notification(node, val, data)
        event_notification(event)

```

hvl_ccb.comm.serial

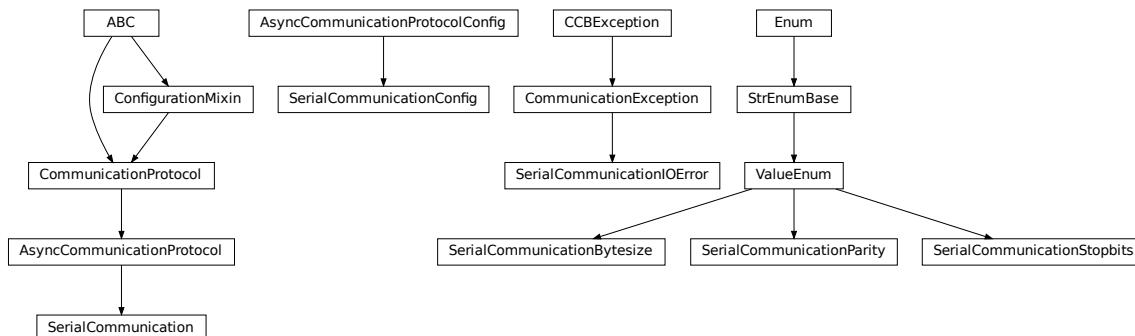

Communication protocol for serial ports. Makes use of the [pySerial](#) library.

```
class SerialCommunication(configuration)
Bases: hvl\_ccb.comm.base.AsyncCommunicationProtocol

Implements the Communication Protocol for serial ports.

close()
    Close the serial connection.

static config_cls()
    Return the default configdataclass class

        Returns a reference to the default configdataclass class

property is_open: bool
    Flag indicating if the serial port is open.

        Returns True if the serial port is open, otherwise False

open()
    Open the serial connection.

        Raises SerialCommunicationIOError – when communication port cannot be opened.

read_bytes() → bytes
    Read the bytes from the serial port till the terminator is found. The input buffer may hold additional lines afterwards.

    This method uses self.access_lock to ensure thread-safety.

        Returns Bytes read from the serial port; b” if there was nothing to read.

        Raises SerialCommunicationIOError – when communication port is not opened

read_single_bytes(size: int = 1) → bytes
    Read the specified number of bytes from the serial port. The input buffer may hold additional data afterwards.

        Returns Bytes read from the serial port; b” if there was nothing to read.

write_bytes(data: bytes) → int
    Write bytes to the serial port.

    This method uses self.access_lock to ensure thread-safety.

        Parameters data – data to write to the serial port

        Returns number of bytes written

        Raises SerialCommunicationIOError – when communication port is not opened

class SerialCommunicationBytesize(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Serial communication bytesize.

EIGHTBITS = 8
FIVEBITS = 5
SEVENBITS = 7
SIXBITS = 6
```

```
class SerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                               encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                               Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,
                               port: Optional[str] = None, baudrate: int = 9600, parity: Union[str,
                               hvl_ccb.comm.serial.SerialCommunicationParity] =
                               SerialCommunicationParity.NONE, stopbits: Union[int, float,
                               hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                               SerialCommunicationStopbits.ONE, bytesize: Union[int,
                               hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                               SerialCommunicationBytesize.EIGHTBITS, timeout: Union[int, float] = 2)
```

Bases: `hvl_ccb.comm.base.AsyncCommunicationProtocolConfig`

Configuration dataclass for `SerialCommunication`.

Bytesize

alias of `hvl_ccb.comm.serial.SerialCommunicationBytesize`

Parity

alias of `hvl_ccb.comm.serial.SerialCommunicationParity`

Stopbits

alias of `hvl_ccb.comm.serial.SerialCommunicationStopbits`

baudrate: int = 9600

Baudrate of the serial port

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

Size of a byte, 5 to 8

clean_values()

create_serial_port() → serial.serialposix.Serial

Create a serial port instance according to specification in this configuration

Returns Closed serial port instance

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

Parity to be used for the connection.

```
port: Optional[str] = None
Port is a string referring to a COM-port (e.g. 'COM3') or a URL. The full list of capabilities is found on the pyserial documentation.
```

```
classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.
```

Returns a list of strings containing all required keys.

```
stopbits: Union[int, float, hvl\_ccb.comm.serial.SerialCommunicationStopbits] = 1
Stopbits setting, can be 1, 1.5 or 2.
```

```
terminator_str() → str
```

```
timeout: Union[int, float] = 2
Timeout in seconds for the serial port
```

```
exception SerialCommunicationIOError
Bases: OSError, hvl\_ccb.comm.base.CommunicationException
Serial communication related I/O errors.
```


```
class SerialCommunicationParity(value=<no_arg>, names=None, module=None, type=None, start=1,
                                  boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum
Serial communication parity.
```

```
EVEN = 'E'
MARK = 'M'
NAMES = {'E': 'Even', 'M': 'Mark', 'N': 'None', 'O': 'Odd', 'S': 'Space'}
NONE = 'N'
ODD = 'O'
SPACE = 'S'
```

```
class SerialCommunicationStopbits(value=<no_arg>, names=None, module=None, type=None, start=1,
                                    boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum
Serial communication stopbits.
```

```
ONE = 1
ONE_POINT_FIVE = 1.5
TWO = 2
```

`hvl_ccb.comm.tcp`

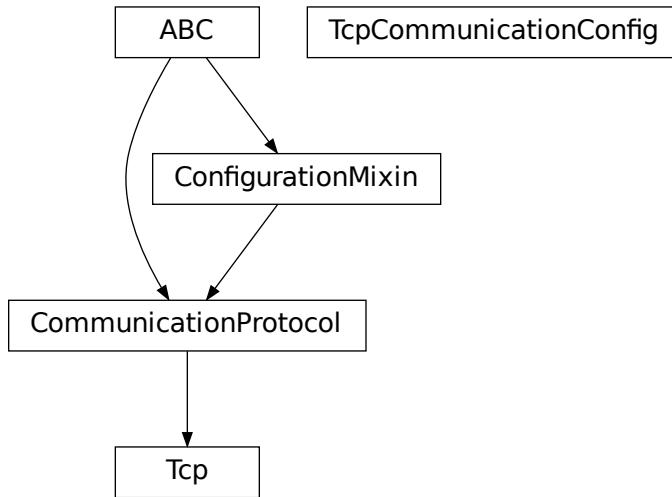

Makes use of the socket library.

TCP communication protocol.

```

class Tcp(configuration)
  Bases: hvl_ccb.comm.base.CommunicationProtocol

  Tcp Communication Protocol.

  close() → None
    Close TCP connection.

  static config_cls() → Type[hvl_ccb.comm.tcp.TcpCommunicationConfig]
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

  open() → None
    Open TCP connection.

  read() → str
    TCP read function :return: information read from TCP buffer formatted as string

  write(command: str = '') → None
    TCP write function :param command: command string to be sent :return: none

class TcpCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], port: int = 54321, bufsize: int = 1024)
  Bases: object

  Configuration dataclass for TcpCommunication.

  bufsize: int = 1024
  clean_values()
  force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.
  
```

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters


- **fieldname** – name of the field
- **value** – value to assign

```
host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]
is_configdataclass = True
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.
    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.
    Returns a list of strings containing all optional keys.

port: int = 54321
classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.
    Returns a list of strings containing all required keys.
```

hvl_ccb.comm.telnet

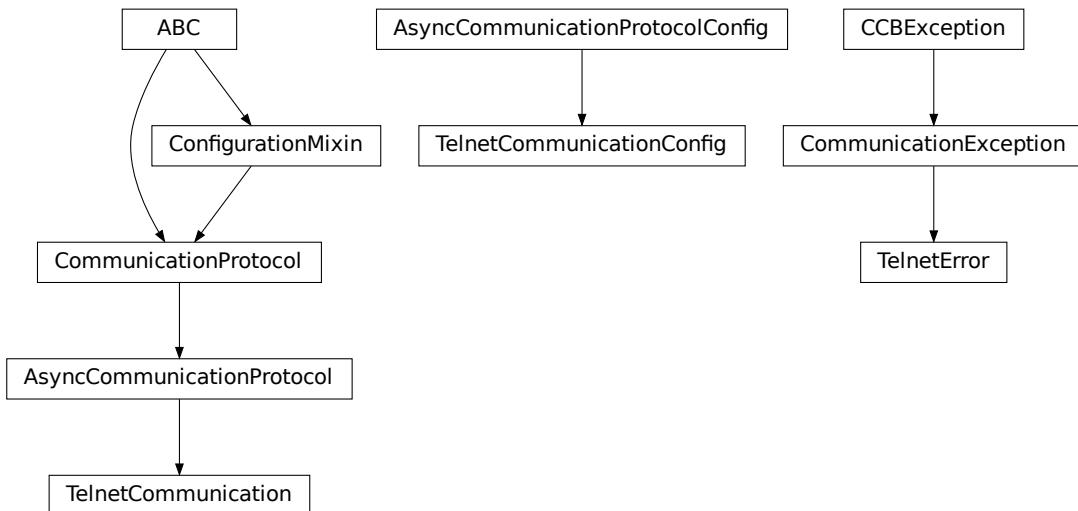

Communication protocol for telnet. Makes use of the `telnetlib` library.

```
class TelnetCommunication(configuration)
Bases: hvl_ccb.comm.base.AsyncCommunicationProtocol
```

Implements the Communication Protocol for telnet.

close()

Close the telnet connection unless it is not closed.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

property is_open: bool

Is the connection open?

Returns True for an open connection

open()

Open the telnet connection unless it is not yet opened.

read_bytes() → bytes

Read data as *bytes* from the telnet connection.

Returns data from telnet connection

Raises *TelnetError* – when connection is not open, raises an Error during the communication

write_bytes(data: bytes)

Write the data as *bytes* to the telnet connection.

Parameters **data** – Data to be sent.

Raises *TelnetError* – when connection is not open, raises an Error during the communication

```
class TelnetCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10,
                                host: Optional[Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]] =
                                None, port: int = 0, timeout: Union[int, float] = 0.2)
```

Bases: *hvl_ccb.comm.base.AsyncCommunicationProtocolConfig*

Configuration dataclass for *TelnetCommunication*.

clean_values()

create_telnet() → Optional[telnetlib.Telnet]

Create a telnet client :return: Opened Telnet object or None if connection is not possible

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: Optional[Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address]] = None

Host to connect to can be localhost or

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 0

Port at which the host is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

timeout: Union[int, float] = 0.2


Timeout for reading a line

exception TelnetError

Bases: OSError, *hvl_ccb.comm.base.CommunicationException*

Telnet communication related errors.

hvl_ccb.comm.visa

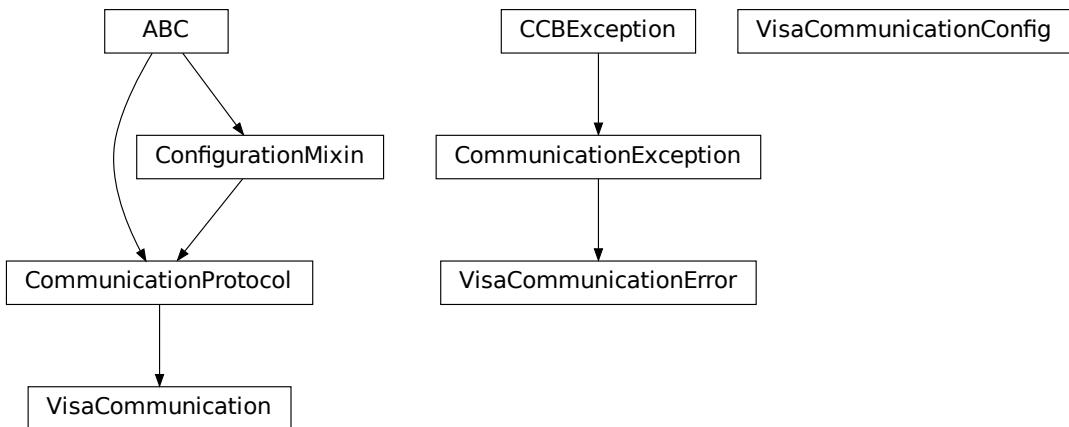

Communication protocol for VISA. Makes use of the pyvisa library. The backend can be NI-Visa or pyvisa-py.

Information on how to install a VISA backend can be found here: https://pyvisa.readthedocs.io/en/master/getting_nivisa.html

So far only TCPIP SOCKET and TCPIP INSTR interfaces are supported.

class VisaCommunication(configuration)

Bases: *hvl_ccb.comm.base.CommunicationProtocol*

Implements the Communication Protocol for VISA / SCPI.

MULTI_COMMANDS_MAX = 5

The maximum of commands that can be sent in one round is 5 according to the VISA standard.

MULTI_COMMANDS_SEPARATOR = ' ; '

The character to separate two commands is ; according to the VISA standard.

WAIT_AFTER_WRITE = 0.08

Small pause in seconds to wait after write operations, allowing devices to really do what we tell them before continuing with further tasks.

close() → None

Close the VISA connection and invalidates the handle.

static config_cls() → Type[*hvl_ccb.comm.visa.VisaCommunicationConfig*]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

open() → None

Open the VISA connection and create the resource.

query(*commands: str) → Union[str, Tuple[str, ...]]

A combination of write(message) and read.

Parameters **commands** – list of commands

Returns list of values

Raises **VisaCommunicationError** – when connection was not started, or when trying to issue too many commands at once.

spoll() → int

Execute serial poll on the device. Reads the status byte register STB. This is a fast function that can be executed periodically in a polling fashion.

Returns integer representation of the status byte

Raises **VisaCommunicationError** – when connection was not started

write(*commands: str) → None

Write commands. No answer is read or expected.

Parameters **commands** – one or more commands to send

Raises **VisaCommunicationError** – when connection was not started

class VisaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], interface_type: Union[str, *hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType*], board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = '')

Bases: object

VISACommunication configuration dataclass.

class InterfaceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: *hvl_ccb.utils.enum.AutoNumberNameEnum*

Supported VISA Interface types.

TCP/IP_INSTR = 2

VXI-11 protocol

TCP/IP_SOCKET = 1

VISA-Raw protocol

address(*host: str, port: Optional[int] = None, board: Optional[int] = None*) → str

Address string specific to the VISA interface type.

Parameters

- **host** – host IP address
- **port** – optional TCP port
- **board** – optional board number

Returns address string

property address: str

Address string depending on the VISA protocol's configuration.

Returns address string corresponding to current configuration

board: int = 0

Board number is typically 0 and comes from old bus systems.

chunk_size: int = 204800

Chunk size is the allocated memory for read operations. The standard is 20kB, and is increased per default here to 200kB. It is specified in bytes.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

host: Union[str, ipAddress.I Pv4Address, ipAddress.I Pv6Address]

interface_type: Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType]

Interface type of the VISA connection, being one of *InterfaceType*.

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

open_timeout: int = 1000

Timeout for opening the connection, in milli seconds.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 5025

TCP port, standard is 5025.

read_termination: str = '\n'

Read termination character.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

timeout: int = 5000

Timeout for commands in milli seconds.

visa_backend: str = ''

Specifies the path to the library to be used with PyVISA as a backend. Defaults to None, which is NI-VISA (if installed), or pyvisa-py (if NI-VISA is not found). To force the use of pyvisa-py, specify '@py' here.

write_termination: str = '\n'

Write termination character.

exception VisaCommunicationError

Bases: OSError, *hvl_ccb.comm.base.CommunicationException*

Base class for VisaCommunication errors.

Module contents

Communication protocols subpackage.


hvl_ccb.dev

Subpackages

hvl_ccb.dev.highland_t560

Submodules

hvl_ccb.dev.highland_t560.base

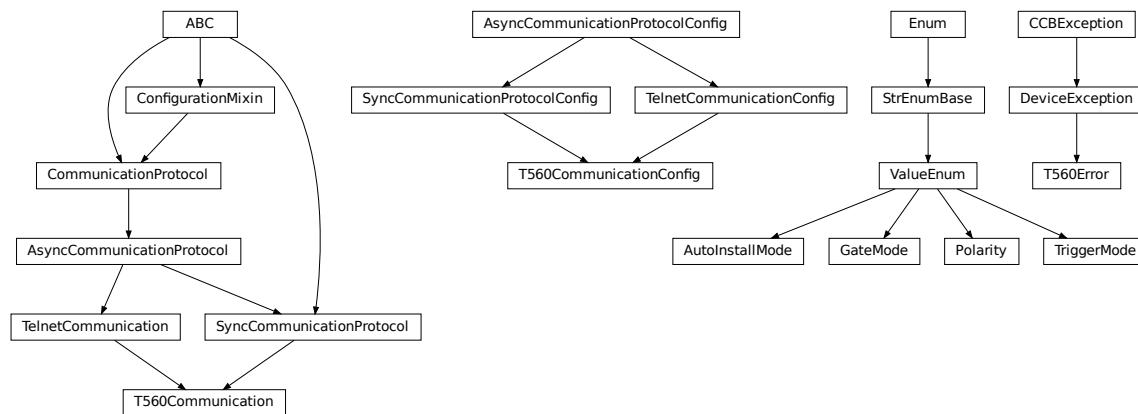

Module containing base device and communication classes and enums.

Communication with device is performed via its ethernet port and a Telnet connection.

class AutoInstallMode(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: *hvl_ccb.utils.enum.ValueEnum*

Modes for installing configuration settings to the device.

INSTALL = 1

OFF = 0

QUEUE = 2

class GateMode(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)
Bases: *hvl_ccb.utils.enum.ValueEnum*

Available T560 gate modes

INPUT = 'INP'

OFF = 'OFF'

OUTPUT = 'OUT'

class Polarity(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)
Bases: *hvl_ccb.utils.enum.ValueEnum*

Possible channel polarity states

ACTIVE_HIGH = 'POS'

ACTIVE_LOW = 'NEG'

class T560Communication(*configuration*)

Bases: *hvl_ccb.comm.base.SyncCommunicationProtocol*, *hvl_ccb.comm.telnet.TelnetCommunication*

Communication class for T560. It uses a TelnetCommunication with the SyncCommunicationProtocol

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(*command: str*) → *str*

Send a command to the device and handle the response.

For device setting queries, response will be 'OK' if successful, or '??' if setting cannot be carried out, raising an error.

Parameters **command** – Command string to be sent

Raises **T560Error** – if no response is received, or if the device responds with an error message.

Returns Response from the device.

class T560CommunicationConfig(*terminator: bytes = b'\r', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address, NoneType] = None, port: int = 2000, timeout: Union[int, float] = 0.2*)

Bases: *hvl_ccb.comm.base.SyncCommunicationProtocolConfig*, *hvl_ccb.comm.telnet.TelnetCommunicationConfig*

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 2000

Port at which the host is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

terminator: bytes = b'\r'

The terminator character. Typically this is b'\r\n' or b'\n', but can also be b'\r' or other combinations.
This defines the end of a single line.

exception T560Error

Bases: *hvl_ccb.dev.base.DeviceException*

T560 related errors.

class TriggerMode(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: *hvl_ccb.utils.enum.ValueEnum*

Available T560 trigger modes

COMMAND = 'REM'

EXT_FALLING_EDGE = 'NEG'


EXT_RISING_EDGE = 'POS'

INT_SYNTHESIZER = 'SYN'

OFF = 'OFF'

hvl_ccb.dev.highland_t560.channel

Module for controlling pulse output channels A, B, C and D.

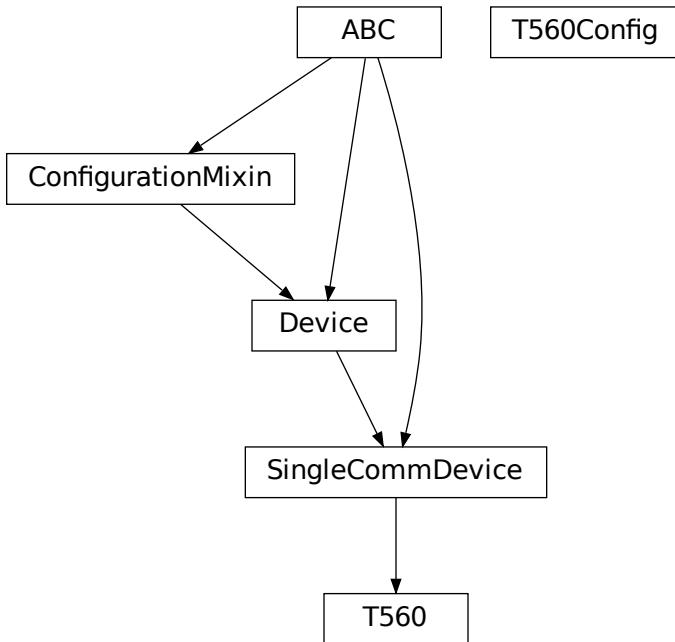hvl_ccb.dev.highland_t560.device

Module for controlling device, including TRIG, CLOCK and GATE I/Os.

```

class T560(com, dev_config=None)
  Bases: hvl_ccb.dev.base.SingleCommDevice

  activate_clock_output()
    Outputs 10 MHz clock signal

  property auto_install_mode: hvl_ccb.dev.highland_t560.base.AutoInstallMode
    Check the autoinstall settings of the T560. The autoinstall mode sets how changes to device settings are applied. See manual section 4.7.2 for more information about these modes.

  property ch_a: hvl_ccb.dev.highland_t560.channel._Channel
    Channel A of T560

  property ch_b: hvl_ccb.dev.highland_t560.channel._Channel
    Channel B of T560

  property ch_c: hvl_ccb.dev.highland_t560.channel._Channel
    Channel C of T560

  property ch_d: hvl_ccb.dev.highland_t560.channel._Channel
    Channel D of T560

  static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class
  
```

```
static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

disarm_trigger()
    Disarm DDG by disabling all trigger sources.

fire_trigger()
    Fire a software trigger.

property frequency: float
    The frequency of the timing cycle in Hz.

property gate_mode: hvl_ccb.dev.highland_t560.base.GateMode
    Check the mode setting of the GATE I/O port.

property gate_polarity: hvl_ccb.dev.highland_t560.base.Polarity
    Check the polarity setting of the GATE I/O port.

load_device_configuration()
    Load the settings saved in nonvolatile memory.

property period: float
    The period of the timing cycle (time between triggers) in seconds.

save_device_configuration()
    Save the current settings to nonvolatile memory.

property trigger_level
    Get external trigger level.

property trigger_mode
    Get device trigger source.

use_external_clock()
    Finds and accepts an external clock signal to the CLOCK input

class T560Config
    Bases: object

    auto_install_mode = 1

    clean_values()
        Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

    force_value(fieldname, value)
        Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.

    Parameters
        • fieldname – name of the field
        • value – value to assign

    is_configdataclass = True

    classmethod keys() → Sequence[str]
        Returns a list of all configdataclass fields key-names.

        Returns a list of strings containing all keys.
```

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

Module contents

This module establishes methods for interfacing with the Highland Technology T560-2 via its ethernet adapter with a telnet communication protocol.

The T560 is a small digital delay & pulse generator. It outputs up to four individually timed pulses with 10-ps precision, given an internal or external trigger.


This module introduces methods for configuring channels, gating, and triggering. Further documentation and a more extensive command list may be obtained from:

<https://www.highlandtechnology.com/DSS/T560DS.shtml>

hvl_ccb.dev.supercube

Submodules

hvl_ccb.dev.supercube.base

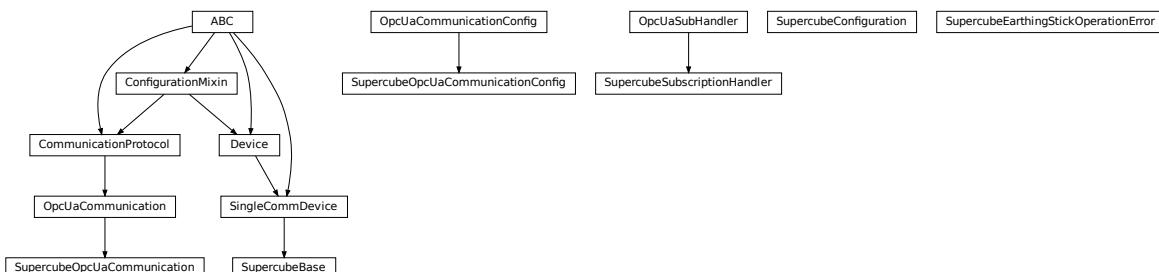

Base classes for the Supercube device.

class SupercubeBase(com, dev_config=None)
Bases: [hvl_ccb.dev.base.SingleCommDevice](#)

Base class for Supercube variants.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

display_message_board() → None

Display 15 newest messages

display_status_board() → None

Display status board.

get_cee16_socket() → bool

Read the on-state of the IEC CEE16 three-phase power socket.

Returns the on-state of the CEE16 power socket

get_door_status(door: int) → *hvl_ccb.dev.supercube.constants.DoorStatus*

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

Parameters `door` – the door number (1..3)

Returns the door status

get_earthing_rod_status(earthing_rod: int) → *hvl_ccb.dev.supercube.constants.EarthRodStatus*

Get the status of a earthing rod. See `constants.EarthRodStatus` for possible returned earthing rod statuses.

Parameters `earthing_rod` – the earthing rod number (1..3)

Returns the earthing rod status

get_earthing_stick_manual(number: int) → *hvl_ccb.dev.supercube.constants.EarthStickOperation*

Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

Parameters `number` – number of the earthing stick (1..6)

Returns operation of the earthing stick in a manual operating mode (open == 0, close == 1)

Raises `ValueError` – when earthing stick number is not valid

get_earthing_stick_operating_status(number: int) →

hvl_ccb.dev.supercube.constants.EarthStickOperatingStatus

Get the operating status of an earthing stick.

Parameters `number` – number of the earthing stick (1..6)

Returns earthing stick operating status (auto == 0, manual == 1)

Raises `ValueError` – when earthing stick number is not valid

get_earthing_stick_status(number: int) → *hvl_ccb.dev.supercube.constants.EarthStickStatus*

Get the status of an earthing stick, whether it is closed, open or undefined (moving).

Parameters `number` – number of the earthing stick (1..6)

Returns earthing stick status

Raises `ValueError` – when earthing stick number is not valid

get_measurement_ratio(channel: int) → float

Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

Parameters `channel` – number of the input channel (1..4)

Returns the ratio

Raises `ValueError` – when channel is not valid

get_measurement_voltage(*channel: int*) → float

Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

Parameters **channel** – number of the input channel (1..4)

Returns measured voltage

Raises **ValueError** – when channel is not valid

get_status() → *hvl_ccb.dev.supercube.constants.SafetyStatus*

Get the safety circuit status of the Supercube. :return: the safety status of the supercube's state machine.

get_support_input(*port: int, contact: int*) → bool

Get the state of a support socket input.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital input read state

Raises **ValueError** – when port or contact number is not valid

get_support_output(*port: int, contact: int*) → bool

Get the state of a support socket output.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital output read state

Raises **ValueError** – when port or contact number is not valid

get_t13_socket(*port: int*) → bool

Read the state of a SEV T13 power socket.

Parameters **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*

Returns on-state of the power socket

Raises **ValueError** – when port is not valid

operate(*state: bool*) → None

Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

Parameters **state** – set operate state

operate_earthing_stick(*number: int, operation:*

hvl_ccb.dev.supercube.constants.EarthingStickOperation) → None

Operation of an earthing stick, which is set to manual operation. If an earthing stick is set to manual, it stays closed even if the system is in states RedReady or RedOperate.

Parameters

- **number** – number of the earthing stick (1..6)
- **operation** – earthing stick manual status (close or open)

Raises **SupercubeEarthingStickOperationError** – when operating status of given number's earthing stick is not manual

quit_error() → None

Quits errors that are active on the Supercube.

read(node_id: str)

Local wrapper for the OPC UA communication protocol read method.

Parameters **node_id** – the id of the node to read.

Returns the value of the variable

ready(state: bool) → None

Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

Parameters **state** – set ready state

set_cee16_socket(state: bool) → None

Switch the IEC CEE16 three-phase power socket on or off.

Parameters **state** – desired on-state of the power socket

Raises **ValueError** – if state is not of type bool

set_message_board(msgs: List[str], display_board: bool = True) → None

Fills messages into message board that display that 15 newest messages with a timestamp.

Parameters

- **msgs** – list of strings
- **display_board** – display 15 newest messages if *True* (default)

Raises **ValueError** – if there are too many messages or the positions indices are invalid.

set_remote_control(state: bool) → None

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

Parameters **state** – desired remote control state

set_status_board(msgs: List[str], pos: Optional[List[int]] = None, clear_board: bool = True, display_board: bool = True) → None

Sets and displays a status board. The messages and the position of the message can be defined.

Parameters

- **msgs** – list of strings
- **pos** – list of integers [0...14]
- **clear_board** – clear unspecified lines if *True* (default), keep otherwise
- **display_board** – display new status board if *True* (default)

Raises **ValueError** – if there are too many messages or the positions indices are invalid.

set_support_output(port: int, contact: int, state: bool) → None

Set the state of a support output socket.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **state** – is the desired state of the support output

Raises **ValueError** – when port or contact number is not valid

set_support_output_impulse(*port: int, contact: int, duration: float = 0.2, pos_pulse: bool = True*) → None

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the *pos_pulse* argument.

This function is blocking.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **duration** – is the length of the impulse in seconds
- **pos_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

Raises ValueError – when port or contact number is not valid

set_t13_socket(*port: int, state: bool*) → None

Set the state of a SEV T13 power socket.

Parameters

- **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*
- **state** – is the desired on-state of the socket

Raises ValueError – when port is not valid or state is not of type bool

start() → None

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

stop() → None

Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

write(*node_id, value*) → None

Local wrapper for the OPC UA communication protocol write method.

Parameters

- **node_id** – the id of the node to read
- **value** – the value to write to the variable

class SupercubeConfiguration(*namespace_index: int = 3, polling_delay_sec: Union[int, float] = 5.0, polling_interval_sec: Union[int, float] = 1.0*)

Bases: object

Configuration dataclass for the Supercube devices.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

```
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

namespace_index: int = 3
    Namespace of the OPC variables, typically this is 3 (coming from Siemens)

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

polling_delay_sec: Union[int, float] = 5.0
polling_interval_sec: Union[int, float] = 1.0

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

exception SupercubeEarthingStickOperationError
    Bases: Exception

class SupercubeOpcUaCommunication(config)
    Bases: hvl_ccb.comm.opc.OpcUaCommunication

    Communication protocol specification for Supercube devices.

    static config_cls()
        Return the default configdataclass class.

        Returns a reference to the default configdataclass class

class SupercubeOpcUaCommunicationConfig(host: Union[str, ipaddress.IPv4Address,
                                                ipaddress.IPv6Address], endpoint_name: str, port: int = 4840,
                                             sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
                                             <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler
                                             object>, update_period: int = 500, wait_timeout_retry_sec:
                                             Union[int, float] = 1, max_timeout_retry_nr: int = 5)
    Bases: hvl_ccb.comm.opc.OpcUaCommunicationConfig

    Communication protocol configuration for OPC UA, specifications for the Supercube devices.

    force_value(fieldname, value)
        Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.

    Parameters
        • fieldname – name of the field
        • value – value to assign

    classmethod keys() → Sequence[str]
        Returns a list of all configdataclass fields key-names.

        Returns a list of strings containing all keys.
```

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

sub_handler: *hvl_ccb.comm.opc.OpcUaSubHandler* =
<*hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler* object>

Subscription handler for data change events

class SupercubeSubscriptionHandler

Bases: *hvl_ccb.comm.opc.OpcUaSubHandler*

OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.


datachange_notification(node: opcua.common.node.Node, val, data)

In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.

Parameters

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

hvl_ccb.dev.supercube.constants

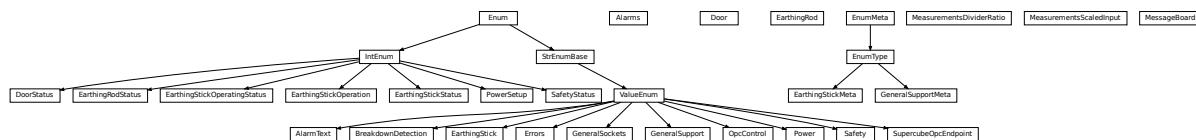

Constants, variable names for the Supercube OPC-connected devices.

class AlarmText(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: *hvl_ccb.utils.enum.ValueEnum*

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the *AlarmText.get()* method to retrieve the enum of an alarm number.

```

Alarm1 = 'STOP Emergency Stop 1'
Alarm10 = 'STOP Earthing stick 2 error while opening'
Alarm11 = 'STOP Earthing stick 3 error while opening'
Alarm12 = 'STOP Earthing stick 4 error while opening'
Alarm13 = 'STOP Earthing stick 5 error while opening'
Alarm14 = 'STOP Earthing stick 6 error while opening'
Alarm15 = 'STOP Earthing stick 1 error while closing'
  
```

```
Alarm16 = 'STOP Earthing stick 2 error while closing'
Alarm17 = 'STOP Earthing stick 3 error while closing'
Alarm18 = 'STOP Earthing stick 4 error while closing'
Alarm19 = 'STOP Earthing stick 5 error while closing'
Alarm2 = 'STOP Emergency Stop 2'
Alarm20 = 'STOP Earthing stick 6 error while closing'
Alarm21 = 'STOP Safety fence 1'
Alarm22 = 'STOP Safety fence 2'
Alarm23 = 'STOP OPC connection error'
Alarm24 = 'STOP Grid power failure'
Alarm25 = 'STOP UPS failure'
Alarm26 = 'STOP 24V PSU failure'
Alarm3 = 'STOP Emergency Stop 3'
Alarm4 = 'STOP Safety Switch 1 error'
Alarm41 = 'WARNING Door 1: Use earthing rod!'
Alarm42 = 'MESSAGE Door 1: Earthing rod is still in setup.'
Alarm43 = 'WARNING Door 2: Use earthing rod!'
Alarm44 = 'MESSAGE Door 2: Earthing rod is still in setup.'
Alarm45 = 'WARNING Door 3: Use earthing rod!'
Alarm46 = 'MESSAGE Door 3: Earthing rod is still in setup.'
Alarm47 = 'MESSAGE UPS charge < 85%'
Alarm48 = 'MESSAGE UPS running on battery'
Alarm5 = 'STOP Safety Switch 2 error'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error while opening'
```

classmethod get(alarm: int)

Get the attribute of this enum for an alarm number.

Parameters `alarm` – the alarm number

Returns the enum for the desired alarm number

not_defined = 'NO ALARM TEXT DEFINED'

```
class Alarms(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._AlarmEnumBase
```

Alarms enumeration containing all variable NodeID strings for the alarm array.

```
Alarm1 = '"DB_Alarm_HMI"."Alarm1"'
```

```
Alarm10 = '"DB_Alarm_HMI"."Alarm10"'
```

```
Alarm100 = '"DB_Alarm_HMI"."Alarm100"'
Alarm101 = '"DB_Alarm_HMI"."Alarm101"'
Alarm102 = '"DB_Alarm_HMI"."Alarm102"'
Alarm103 = '"DB_Alarm_HMI"."Alarm103"'
Alarm104 = '"DB_Alarm_HMI"."Alarm104"'
Alarm105 = '"DB_Alarm_HMI"."Alarm105"'
Alarm106 = '"DB_Alarm_HMI"."Alarm106"'
Alarm107 = '"DB_Alarm_HMI"."Alarm107"'
Alarm108 = '"DB_Alarm_HMI"."Alarm108"'
Alarm109 = '"DB_Alarm_HMI"."Alarm109"'
Alarm11 = '"DB_Alarm_HMI"."Alarm11"'
Alarm110 = '"DB_Alarm_HMI"."Alarm110"'
Alarm111 = '"DB_Alarm_HMI"."Alarm111"'
Alarm112 = '"DB_Alarm_HMI"."Alarm112"'
Alarm113 = '"DB_Alarm_HMI"."Alarm113"'
Alarm114 = '"DB_Alarm_HMI"."Alarm114"'
Alarm115 = '"DB_Alarm_HMI"."Alarm115"'
Alarm116 = '"DB_Alarm_HMI"."Alarm116"'
Alarm117 = '"DB_Alarm_HMI"."Alarm117"'
Alarm118 = '"DB_Alarm_HMI"."Alarm118"'
Alarm119 = '"DB_Alarm_HMI"."Alarm119"'
Alarm12 = '"DB_Alarm_HMI"."Alarm12"'
Alarm120 = '"DB_Alarm_HMI"."Alarm120"'
Alarm121 = '"DB_Alarm_HMI"."Alarm121"'
Alarm122 = '"DB_Alarm_HMI"."Alarm122"'
Alarm123 = '"DB_Alarm_HMI"."Alarm123"'
Alarm124 = '"DB_Alarm_HMI"."Alarm124"'
Alarm125 = '"DB_Alarm_HMI"."Alarm125"'
Alarm126 = '"DB_Alarm_HMI"."Alarm126"'
Alarm127 = '"DB_Alarm_HMI"."Alarm127"'
Alarm128 = '"DB_Alarm_HMI"."Alarm128"'
Alarm129 = '"DB_Alarm_HMI"."Alarm129"'
Alarm13 = '"DB_Alarm_HMI"."Alarm13"'
Alarm130 = '"DB_Alarm_HMI"."Alarm130"'
Alarm131 = '"DB_Alarm_HMI"."Alarm131"'
Alarm132 = '"DB_Alarm_HMI"."Alarm132"'
```

```
Alarm133 = '"DB_Alarm_HMI"."Alarm133"'
Alarm134 = '"DB_Alarm_HMI"."Alarm134"'
Alarm135 = '"DB_Alarm_HMI"."Alarm135"'
Alarm136 = '"DB_Alarm_HMI"."Alarm136"'
Alarm137 = '"DB_Alarm_HMI"."Alarm137"'
Alarm138 = '"DB_Alarm_HMI"."Alarm138"'
Alarm139 = '"DB_Alarm_HMI"."Alarm139"'
Alarm14 = '"DB_Alarm_HMI"."Alarm14"'
Alarm140 = '"DB_Alarm_HMI"."Alarm140"'
Alarm141 = '"DB_Alarm_HMI"."Alarm141"'
Alarm142 = '"DB_Alarm_HMI"."Alarm142"'
Alarm143 = '"DB_Alarm_HMI"."Alarm143"'
Alarm144 = '"DB_Alarm_HMI"."Alarm144"'
Alarm145 = '"DB_Alarm_HMI"."Alarm145"'
Alarm146 = '"DB_Alarm_HMI"."Alarm146"'
Alarm147 = '"DB_Alarm_HMI"."Alarm147"'
Alarm148 = '"DB_Alarm_HMI"."Alarm148"'
Alarm149 = '"DB_Alarm_HMI"."Alarm149"'
Alarm15 = '"DB_Alarm_HMI"."Alarm15"'
Alarm150 = '"DB_Alarm_HMI"."Alarm150"'
Alarm151 = '"DB_Alarm_HMI"."Alarm151"'
Alarm16 = '"DB_Alarm_HMI"."Alarm16"'
Alarm17 = '"DB_Alarm_HMI"."Alarm17"'
Alarm18 = '"DB_Alarm_HMI"."Alarm18"'
Alarm19 = '"DB_Alarm_HMI"."Alarm19"'
Alarm2 = '"DB_Alarm_HMI"."Alarm2"'
Alarm20 = '"DB_Alarm_HMI"."Alarm20"'
Alarm21 = '"DB_Alarm_HMI"."Alarm21"'
Alarm22 = '"DB_Alarm_HMI"."Alarm22"'
Alarm23 = '"DB_Alarm_HMI"."Alarm23"'
Alarm24 = '"DB_Alarm_HMI"."Alarm24"'
Alarm25 = '"DB_Alarm_HMI"."Alarm25"'
Alarm26 = '"DB_Alarm_HMI"."Alarm26"'
Alarm27 = '"DB_Alarm_HMI"."Alarm27"'
Alarm28 = '"DB_Alarm_HMI"."Alarm28"'
Alarm29 = '"DB_Alarm_HMI"."Alarm29"'
```

```
Alarm3 = '"DB_Alarm_HMI"."Alarm3"'
Alarm30 = '"DB_Alarm_HMI"."Alarm30"'
Alarm31 = '"DB_Alarm_HMI"."Alarm31"'
Alarm32 = '"DB_Alarm_HMI"."Alarm32"'
Alarm33 = '"DB_Alarm_HMI"."Alarm33"'
Alarm34 = '"DB_Alarm_HMI"."Alarm34"'
Alarm35 = '"DB_Alarm_HMI"."Alarm35"'
Alarm36 = '"DB_Alarm_HMI"."Alarm36"'
Alarm37 = '"DB_Alarm_HMI"."Alarm37"'
Alarm38 = '"DB_Alarm_HMI"."Alarm38"'
Alarm39 = '"DB_Alarm_HMI"."Alarm39"'
Alarm4 = '"DB_Alarm_HMI"."Alarm4"'
Alarm40 = '"DB_Alarm_HMI"."Alarm40"'
Alarm41 = '"DB_Alarm_HMI"."Alarm41"'
Alarm42 = '"DB_Alarm_HMI"."Alarm42"'
Alarm43 = '"DB_Alarm_HMI"."Alarm43"'
Alarm44 = '"DB_Alarm_HMI"."Alarm44"'
Alarm45 = '"DB_Alarm_HMI"."Alarm45"'
Alarm46 = '"DB_Alarm_HMI"."Alarm46"'
Alarm47 = '"DB_Alarm_HMI"."Alarm47"'
Alarm48 = '"DB_Alarm_HMI"."Alarm48"'
Alarm49 = '"DB_Alarm_HMI"."Alarm49"'
Alarm5 = '"DB_Alarm_HMI"."Alarm5"'
Alarm50 = '"DB_Alarm_HMI"."Alarm50"'
Alarm51 = '"DB_Alarm_HMI"."Alarm51"'
Alarm52 = '"DB_Alarm_HMI"."Alarm52"'
Alarm53 = '"DB_Alarm_HMI"."Alarm53"'
Alarm54 = '"DB_Alarm_HMI"."Alarm54"'
Alarm55 = '"DB_Alarm_HMI"."Alarm55"'
Alarm56 = '"DB_Alarm_HMI"."Alarm56"'
Alarm57 = '"DB_Alarm_HMI"."Alarm57"'
Alarm58 = '"DB_Alarm_HMI"."Alarm58"'
Alarm59 = '"DB_Alarm_HMI"."Alarm59"'
Alarm6 = '"DB_Alarm_HMI"."Alarm6"'
Alarm60 = '"DB_Alarm_HMI"."Alarm60"'
Alarm61 = '"DB_Alarm_HMI"."Alarm61"'
```

```
Alarm62 = '"DB_Alarm_HMI"."Alarm62"'
Alarm63 = '"DB_Alarm_HMI"."Alarm63"'
Alarm64 = '"DB_Alarm_HMI"."Alarm64"'
Alarm65 = '"DB_Alarm_HMI"."Alarm65"'
Alarm66 = '"DB_Alarm_HMI"."Alarm66"'
Alarm67 = '"DB_Alarm_HMI"."Alarm67"'
Alarm68 = '"DB_Alarm_HMI"."Alarm68"'
Alarm69 = '"DB_Alarm_HMI"."Alarm69"'
Alarm7 = '"DB_Alarm_HMI"."Alarm7"'
Alarm70 = '"DB_Alarm_HMI"."Alarm70"'
Alarm71 = '"DB_Alarm_HMI"."Alarm71"'
Alarm72 = '"DB_Alarm_HMI"."Alarm72"'
Alarm73 = '"DB_Alarm_HMI"."Alarm73"'
Alarm74 = '"DB_Alarm_HMI"."Alarm74"'
Alarm75 = '"DB_Alarm_HMI"."Alarm75"'
Alarm76 = '"DB_Alarm_HMI"."Alarm76"'
Alarm77 = '"DB_Alarm_HMI"."Alarm77"'
Alarm78 = '"DB_Alarm_HMI"."Alarm78"'
Alarm79 = '"DB_Alarm_HMI"."Alarm79"'
Alarm8 = '"DB_Alarm_HMI"."Alarm8"'
Alarm80 = '"DB_Alarm_HMI"."Alarm80"'
Alarm81 = '"DB_Alarm_HMI"."Alarm81"'
Alarm82 = '"DB_Alarm_HMI"."Alarm82"'
Alarm83 = '"DB_Alarm_HMI"."Alarm83"'
Alarm84 = '"DB_Alarm_HMI"."Alarm84"'
Alarm85 = '"DB_Alarm_HMI"."Alarm85"'
Alarm86 = '"DB_Alarm_HMI"."Alarm86"'
Alarm87 = '"DB_Alarm_HMI"."Alarm87"'
Alarm88 = '"DB_Alarm_HMI"."Alarm88"'
Alarm89 = '"DB_Alarm_HMI"."Alarm89"'
Alarm9 = '"DB_Alarm_HMI"."Alarm9"'
Alarm90 = '"DB_Alarm_HMI"."Alarm90"'
Alarm91 = '"DB_Alarm_HMI"."Alarm91"'
Alarm92 = '"DB_Alarm_HMI"."Alarm92"'
Alarm93 = '"DB_Alarm_HMI"."Alarm93"'
Alarm94 = '"DB_Alarm_HMI"."Alarm94"'
```

```

Alarm95 = '"DB_Alarm_HMI"."Alarm95"'
Alarm96 = '"DB_Alarm_HMI"."Alarm96"'
Alarm97 = '"DB_Alarm_HMI"."Alarm97"'
Alarm98 = '"DB_Alarm_HMI"."Alarm98"'
Alarm99 = '"DB_Alarm_HMI"."Alarm99"'

class BreakdownDetection(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Node ID strings for the breakdown detection.

TODO: these variable NodeIDs are not tested and/or correct yet.

activated = '"Ix_Allg_Breakdown_activated"'
Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the system or not.

reset = '"Qx_Allg_Breakdown_reset"'
Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

triggered = '"Ix_Allg_Breakdown_triggered"'
Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using the safety circuit state, therefore no method is implemented to read this out directly.

class Door(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._DoorEnumBase

Variable NodeID strings for doors.

status_1 = '"DB_Safety_Circuit"."Door_1"."si_HMI_status"'
status_2 = '"DB_Safety_Circuit"."Door_2"."si_HMI_status"'
status_3 = '"DB_Safety_Circuit"."Door_3"."si_HMI_status"'

class DoorStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for doors.

closed = 2
Door is closed, but not locked.

error = 4
Door has an error or was opened in locked state (either with emergency stop or from the inside).

inactive = 0
not enabled in Supercube HMI setup, this door is not supervised.

locked = 3
Door is closed and locked (safe state).

open = 1
Door is open.

class EarthingRod(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.dev.supercube.constants._DoorEnumBase

Variable NodeID strings for earthing rods.

status_1 = '"DB_Safety_Circuit"."Door_1"."Ix_earthingrod"'

```

```
status_2 = '"DB_Safety_Circuit"."Door_2"."Ix_earthingrod"'
status_3 = '"DB_Safety_Circuit"."Door_3"."Ix_earthingrod"'

class EarthingRodStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
                        boundary=None)
Bases: aenum.IntEnum

Possible status values for earthing rods.

experiment_blocked = 0
    earthing rod is somewhere in the experiment and blocks the start of the experiment

experiment_ready = 1
    earthing rod is hanging next to the door, experiment is ready to operate

class EarthingStick(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the
earthing in manual mode.

classmethod manual(number: int)
    Get the manual enum instance for an earthing stick number.

    Parameters number – the earthing stick (1..6)

    Returns the manual instance

    Raises ValueError – when earthing stick number is not valid

manual_1 = '"DB_Safety_Circuit"."Earthstick_1"."sx_earthing_manually"'
manual_2 = '"DB_Safety_Circuit"."Earthstick_2"."sx_earthing_manually"'
manual_3 = '"DB_Safety_Circuit"."Earthstick_3"."sx_earthing_manually"'
manual_4 = '"DB_Safety_Circuit"."Earthstick_4"."sx_earthing_manually"'
manual_5 = '"DB_Safety_Circuit"."Earthstick_5"."sx_earthing_manually"'
manual_6 = '"DB_Safety_Circuit"."Earthstick_6"."sx_earthing_manually"'

classmethod manuals() → Tuple[hvl_ccb.dev.supercube.constants.EarthlingStick, ...]
    Get all earthing stick manual instances.

    Returns tuple of manual instances

property number: int
    Get corresponding earthing stick number.

    Returns earthing stick number (1..6)

classmethod operating_status(number: int)
    Get the operating status enum instance for an earthing stick number.

    Parameters number – the earthing stick (1..6)

    Returns the operating status instance

    Raises ValueError – when earthing stick number is not valid

operating_status_1 = '"DB_Safety_Circuit"."Earthstick_1"."sx_manual_control_active"'
operating_status_2 = '"DB_Safety_Circuit"."Earthstick_2"."sx_manual_control_active"'
operating_status_3 = '"DB_Safety_Circuit"."Earthstick_3"."sx_manual_control_active"'
```

```
operating_status_4 = '"DB_Safety_Circuit"."Earthstick_4"."sx_manual_control_active"'
operating_status_5 = '"DB_Safety_Circuit"."Earthstick_5"."sx_manual_control_active"'
operating_status_6 = '"DB_Safety_Circuit"."Earthstick_6"."sx_manual_control_active"'
classmethod operating_statuses() → Tuple[hvl_ccb.dev.supercube.constants.EarthingStick, ...]
```

Get all earthing stick operating status instances.

Returns tuple of operating status instances

```
classmethod range() → Sequence[int]
```

Integer range of all earthing sticks.

Returns sequence of earthing sticks numbers

```
classmethod status(number: int) → hvl_ccb.dev.supercube.constants.EarthingStick
```

Get the status enum instance for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status instance

Raises **ValueError** – when earthing stick number is not valid

```
status_1 = '"DB_Safety_Circuit"."Earthstick_1"."si_HMI_Status"'
status_2 = '"DB_Safety_Circuit"."Earthstick_2"."si_HMI_Status"'
status_3 = '"DB_Safety_Circuit"."Earthstick_3"."si_HMI_Status"'
status_4 = '"DB_Safety_Circuit"."Earthstick_4"."si_HMI_Status"'
status_5 = '"DB_Safety_Circuit"."Earthstick_5"."si_HMI_Status"'
status_6 = '"DB_Safety_Circuit"."Earthstick_6"."si_HMI_Status"
```

```
classmethod statuses() → Tuple[hvl_ccb.dev.supercube.constants.EarthingStick, ...]
```

Get all earthing stick status instances.

Returns tuple of status instances

```
class EarthingStickMeta(clsname, bases, clsdic, **kwargs)
```

Bases: aenum.EnumType

```
class EarthingStickOperatingStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
                                    boundary=None)
```

Bases: aenum.IntEnum

Operating Status for an earthing stick. Stick can be used in auto or manual mode.

auto = 0

manual = 1

```
class EarthingStickOperation(value=<no_arg>, names=None, module=None, type=None, start=1,
                                boundary=None)
```

Bases: aenum.IntEnum

Operation of the earthing stick in manual operating mode. Can be closed or opened.

close = 1

open = 0

```
class EarthingStickStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
Bases: aenum.IntEnum
Status of an earthing stick. These are the possible values in the status integer e.g. in EarthingStick.status_1.
```

closed = 1
Earthing is closed (safe).

error = 3
Earthing is in error, e.g. when the stick did not close correctly or could not open.

inactive = 0
Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be enabled in the Supercube HMI setup.

open = 2
Earthing is open (not safe).

```
class Errors(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
Variable NodeID strings for information regarding error, warning and message handling.
```

message = '"DB_Message_Buffer"."Info_active"
Boolean read-only variable telling if a message is active.

quit = '"DB_Message_Buffer"."Reset_button"
Writable boolean for the error quit button.

stop = '"DB_Message_Buffer"."Stop_active"
Boolean read-only variable telling if a stop is active.

warning = '"DB_Message_Buffer"."Warning_active"
Boolean read-only variable telling if a warning is active.

```
class GeneralSockets(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
NodeID strings for the power sockets (3x T13 and 1xCEE16).
```

cee16 = '"Qx_Allg_Socket_CEE16"
CEE16 socket (writeable boolean).

t13_1 = '"Qx_Allg_Socket_T13_1"
SEV T13 socket No. 1 (writable boolean).

t13_2 = '"Qx_Allg_Socket_T13_2"
SEV T13 socket No. 2 (writable boolean).

t13_3 = '"Qx_Allg_Socket_T13_3"
SEV T13 socket No. 3 (writable boolean).

```
class GeneralSupport(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
NodeID strings for the support inputs and outputs.
```

classmethod contact_range() → Sequence[int]
Integer range of all contacts.

Returns sequence of contact numbers

in_1_1 = '"Ix_Allg_Support1_1"

```

in_1_2 = '"Ix_Allg_Support1_2"'
in_2_1 = '"Ix_Allg_Support2_1"'
in_2_2 = '"Ix_Allg_Support2_2"'
in_3_1 = '"Ix_Allg_Support3_1"'
in_3_2 = '"Ix_Allg_Support3_2"'
in_4_1 = '"Ix_Allg_Support4_1"'
in_4_2 = '"Ix_Allg_Support4_2"'
in_5_1 = '"Ix_Allg_Support5_1"'
in_5_2 = '"Ix_Allg_Support5_2"'
in_6_1 = '"Ix_Allg_Support6_1"'
in_6_2 = '"Ix_Allg_Support6_2"'

classmethod input(port: int, contact: int)
    Get the NodeID string for a support input.

```

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string**Raises ValueError** – when port or contact number is not valid

```

out_1_1 = '"Qx_Allg_Support1_1"'
out_1_2 = '"Qx_Allg_Support1_2"'
out_2_1 = '"Qx_Allg_Support2_1"'
out_2_2 = '"Qx_Allg_Support2_2"'
out_3_1 = '"Qx_Allg_Support3_1"'
out_3_2 = '"Qx_Allg_Support3_2"'
out_4_1 = '"Qx_Allg_Support4_1"'
out_4_2 = '"Qx_Allg_Support4_2"'
out_5_1 = '"Qx_Allg_Support5_1"'
out_5_2 = '"Qx_Allg_Support5_2"'
out_6_1 = '"Qx_Allg_Support6_1"'
out_6_2 = '"Qx_Allg_Support6_2"'

classmethod output(port: int, contact: int)
    Get the NodeID string for a support output.

```

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string**Raises ValueError** – when port or contact number is not valid

```
classmethod port_range() → Sequence[int]
```

Integer range of all ports.

Returns sequence of port numbers

```
class GeneralSupportMeta(clsname, bases, clsdic, **kwargs)
```

Bases: aenum.EnumType

```
class MeasurementsDividerRatio(value=<no_arg>, names=None, module=None, type=None, start=1,  
                                boundary=None)
```

Bases: hvl_ccb.dev.supercube.constants._InputEnumBase

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

```
input_1 = '"DB_Measurements"."si_Divider_Ratio_1"'
```

```
input_2 = '"DB_Measurements"."si_Divider_Ratio_2"'
```

```
input_3 = '"DB_Measurements"."si_Divider_Ratio_3"'
```

```
input_4 = '"DB_Measurements"."si_Divider_Ratio_4"'
```

```
class MeasurementsScaledInput(value=<no_arg>, names=None, module=None, type=None, start=1,  
                                boundary=None)
```

Bases: hvl_ccb.dev.supercube.constants._InputEnumBase

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in [MeasurementsDividerRatio](#).

```
input_1 = '"DB_Measurements"."si_scaled_Voltage_Input_1"'
```

```
input_2 = '"DB_Measurements"."si_scaled_Voltage_Input_2"'
```

```
input_3 = '"DB_Measurements"."si_scaled_Voltage_Input_3"'
```

```
input_4 = '"DB_Measurements"."si_scaled_Voltage_Input_4"'
```

```
class MessageBoard(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: hvl_ccb.dev.supercube.constants._LineEnumBase

Variable NodeID strings for message board lines.

```
line_1 = '"DB_OP_C_Connection"."Is_status_Line_1"'
```

```
line_10 = '"DB_OP_C_Connection"."Is_status_Line_10"'
```

```
line_11 = '"DB_OP_C_Connection"."Is_status_Line_11"'
```

```
line_12 = '"DB_OP_C_Connection"."Is_status_Line_12"'
```

```
line_13 = '"DB_OP_C_Connection"."Is_status_Line_13"'
```

```
line_14 = '"DB_OP_C_Connection"."Is_status_Line_14"'
```

```
line_15 = '"DB_OP_C_Connection"."Is_status_Line_15"'
```

```
line_2 = '"DB_OP_C_Connection"."Is_status_Line_2"'
```

```
line_3 = '"DB_OP_C_Connection"."Is_status_Line_3"'
```

```
line_4 = '"DB_OP_C_Connection"."Is_status_Line_4"'
```

```
line_5 = '"DB_OP_C_Connection"."Is_status_Line_5"'
```

```
line_6 = '"DB_OP_C_Connection"."Is_status_Line_6"'
```

```

line_7 = '"DB_OPCT_Connection"."Is_status_Line_7"'
line_8 = '"DB_OPCT_Connection"."Is_status_Line_8"'
line_9 = '"DB_OPCT_Connection"."Is_status_Line_9"'

class OpcControl(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Variable NodeID strings for supervision of the OPC connection from the controlling workstation to the Supercube.

active = '"DB_OPCT_Connection"."sx_OPCT_active"'
writable boolean to enable OPC remote control and display a message window on the Supercube HMI.

live = '"DB_OPCT_Connection"."sx_OPCT_lifebit"'

class Power(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl\_ccb.utils.enum.ValueEnum

Variable NodeID strings concerning power data.

TODO: these variable NodeIDs are not tested and/or correct yet, they don't exist yet on Supercube side.

current_primary = 'Qr_Power_FU_actual_Current'
Primary current in ampere, measured by the frequency converter. (read-only)

frequency = 'Ir_Power_FU_Frequency'
Frequency converter output frequency. (read-only)

setup = 'Qi_Power_Setup'
Power setup that is configured using the Supercube HMI. The value corresponds to the ones in PowerSetup. (read-only)

voltage_max = 'Iw_Power_max_Voltage'
Maximum voltage allowed by the current experimental setup. (read-only)

voltage_primary = 'Qr_Power_FU_actual_Voltage'
Primary voltage in volts, measured by the frequency converter at its output. (read-only)

voltage_slope = 'Ir_Power_dUDt'
Voltage slope in V/s.

voltage_target = 'Ir_Power_Target_Voltage'
Target voltage setpoint in V.

class PowerSetup(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible power setups corresponding to the value of variable Power.setup.

AC_DoubleStage_150kV = 4
AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum voltage of 150kV.

AC_DoubleStage_200kV = 5
AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

AC_SingleStage_100kV = 3
AC voltage with MWB transformer set to 100kV maximum voltage.

AC_SingleStage_50kV = 2
AC voltage with MWB transformer set to 50kV maximum voltage.

```

DC_DoubleStage_280kV = 8

DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single stage transformer with Greinacher voltage doubling rectifier)

DC_SingleStage_140kV = 7

DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

External = 1

External power supply fed through blue CEE32 input using isolation transformer and safety switches of the Supercube, or using an external safety switch attached to the Supercube Type B.

Internal = 6

Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

NoPower = 0

No safety switches, use only safety components (doors, fence, earthing...) without any power.

class Safety(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: `hvl_ccb.utils.enum.ValueEnum`

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

status = "DB_Safety_Circuit"."si_safe_status"

Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in `SafetyStatus`.

switch_to_operate = "DB_Safety_Circuit"."sx_safe_switch_to_operate"

Writable boolean for switching to Red Operate (locket, HV on) state.

switch_to_ready = "DB_Safety_Circuit"."sx_safe_switch_to_ready"

Writable boolean for switching to Red Ready (locked, HV off) state.

class SafetyStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: `aenum.IntEnum`

Safety status values that are possible states returned from `hvl_ccb.dev.supercube.base.Supercube.get_status()`. These values correspond to the states of the Supercube’s safety circuit statemachine.

Error = 6

System is in error mode.

GreenNotReady = 1

System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

GreenReady = 2

System is safe and all safety elements are in place to be able to switch to *ready*.

Initializing = 0

System is initializing or booting.

QuickStop = 5

Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

RedOperate = 4

System is locked in red state and in *operate* mode, i.e. high voltage on.

RedReady = 3

System is locked in red state and *ready* to go to *operate* mode.

```
class SupercubeOpcEndpoint(value=<no_arg>, names=None, module=None, type=None, start=1,
                             boundary=None)
```

Bases: `hvl_ccb.utils.enum.ValueEnum`

OPC Server Endpoint strings for the supercube variants.


`A = 'Supercube Typ A'`

`B = 'Supercube Typ B'`

`T13_SOCKET_PORTS = (1, 2, 3)`

Port numbers of SEV T13 power socket

`hvl_ccb.dev.supercube.typ_a`

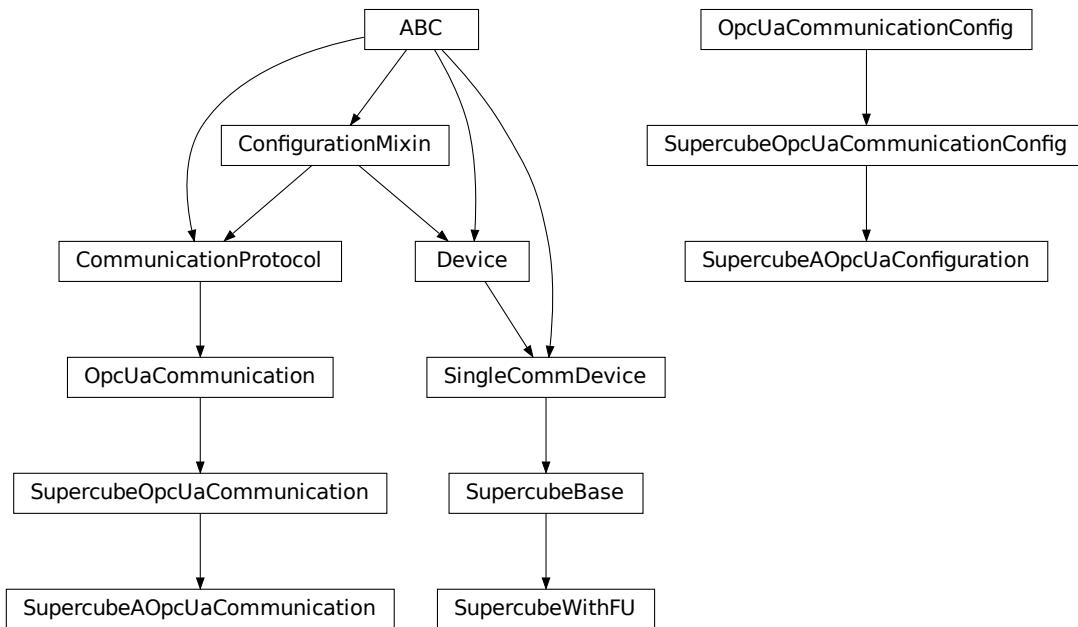

Supercube Typ A module.

```
class SupercubeAOpcUaCommunication(config)
```

Bases: `hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunication`

`static config_cls()`

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class SupercubeAOpcUaConfiguration(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address],
                                    endpoint_name: str = 'Supercube Typ A', port: int = 4840,
                                    sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =
                                    <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object at
                                    0x7fb04aa95850>, update_period: int = 500, wait_timeout_retry_sec:
                                    Union[int, float] = 1, max_timeout_retry_nr: int = 5)
```

Bases: `hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunicationConfig`

```
endpoint_name: str = 'Supercube Typ A'  
Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'  
force_value(fieldname, value)  
Forces a value to a dataclass field despite the class being frozen.  
NOTE: you can define post_force_value method with same signature as this method to do extra processing  
after value has been forced on fieldname.  
Parameters

- fieldname – name of the field
- value – value to assign

classmethod keys() → Sequence[str]  
Returns a list of all configdataclass fields key-names.  
Returns a list of strings containing all keys.  
classmethod optional_defaults() → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified  
on instantiation.  
Returns a list of strings containing all optional keys.  
classmethod required_keys() → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on  
instantiation.  
Returns a list of strings containing all required keys.  
class SupercubeWithFU(com, dev_config=None)  
Bases: hvl\_ccb.dev.supercube.base.SuperCubeBase  
Variant A of the Supercube with frequency converter.  
static default_com_cls()  
Get the class for the default communication protocol used with this device.  
Returns the type of the standard communication protocol for this device  
fso_reset() → None  
TODO: test fso_reset with device  
Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.  
get_frequency() → float  
TODO: test get_frequency with device  
Read the electrical frequency of the current Supercube setup.  
Returns the frequency in Hz  
get_fso_active() → bool  
TODO: test get_fso_active with device  
Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.  
Returns state of the FSO functionality  
get_max_voltage() → float  
TODO: test get_max_voltage with device  
Reads the maximum voltage of the setup and returns in V.  
Returns the maximum voltage of the setup in V.
```

get_power_setup() → *hvl_ccb.dev.supercube.constants.PowerSetup*

TODO: test **get_power_setup** with device

Return the power setup selected in the Supercube's settings.

Returns the power setup

get_primary_current() → float

TODO: test **get_primary_current** with device

Read the current primary current at the output of the frequency converter (before transformer).

Returns primary current in A

get_primary_voltage() → float

TODO: test **get_primary_voltage** with device

Read the current primary voltage at the output of the frequency converter (before transformer).

Returns primary voltage in V

get_target_voltage() → float

TODO: test **get_target_voltage** with device

Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to **set_target_voltage()**.

Returns the setpoint voltage in V.

set_slope(*slope: float*) → None

TODO: test **set_slope** with device

Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.

Parameters **slope** – voltage slope in V/s (0..15)


set_target_voltage(*volt_v: float*) → None

TODO: test **set_target_voltage** with device

Set the output voltage to a defined value in V.

Parameters **volt_v** – the desired voltage in V

`hvl_ccb.dev.supercube.typ_b`

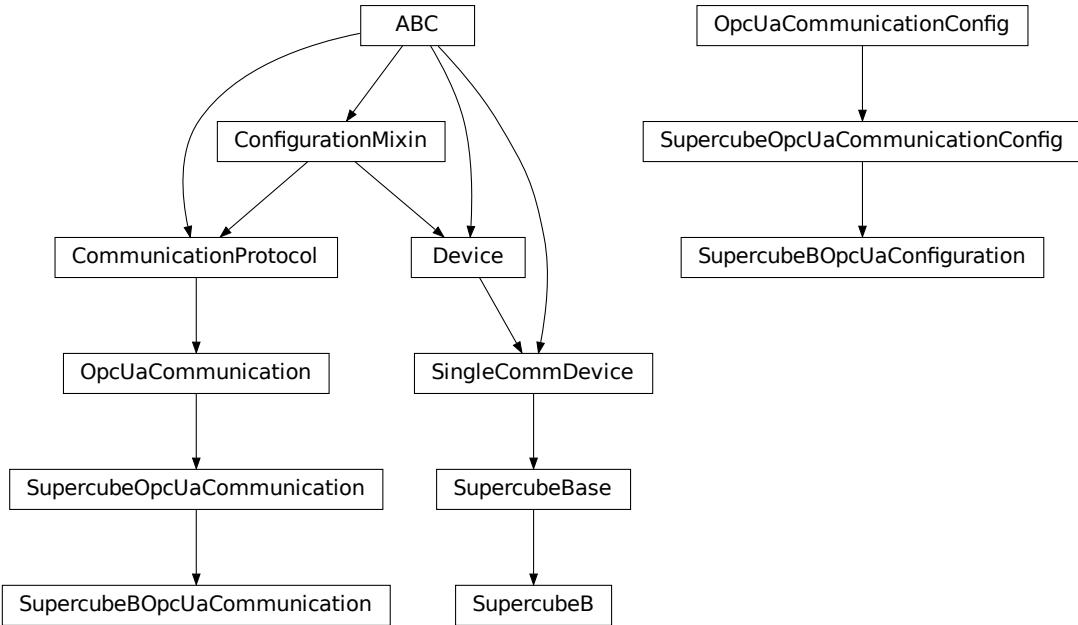

Supercube Typ B module.

class SupercubeB(*com, dev_config=None*)
Bases: *hvl_ccb.dev.supercube.base.SupercubeBase*

Variant B of the Supercube without frequency converter but external safety switches.

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

class SupercubeBOpcaCommunication(*config*)
Bases: *hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunication*

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class SupercubeBOpcaConfiguration(*host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], endpoint_name: str = 'Supercube Typ B', port: int = 4840, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube.base.SupercubeSubscriptionHandler object at 0x7fb04aa95850>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5*)
Bases: *hvl_ccb.dev.supercube.base.SupercubeOpcUaCommunicationConfig*

endpoint_name: str = 'Supercube Typ B'

Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.


classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

Module contents

Supercube package with implementation for system versions from 2019 on (new concept with hard-PLC Siemens S7-1500 as CPU).

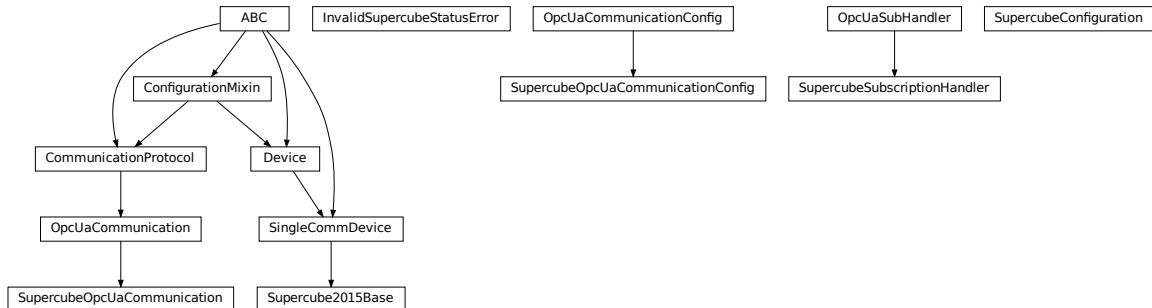hvl_ccb.dev.supercube2015**Submodules****hvl_ccb.dev.supercube2015.base**

Base classes for the Supercube device.

exception InvalidSupercubeStatusError

Bases: `Exception`

Exception raised when supercube has invalid status.

class Supercube2015Base(*com, dev_config=None*)

Bases: `hvl_ccb.dev.base.SingleCommDevice`

Base class for Supercube variants.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_cee16_socket() → bool

Read the on-state of the IEC CEE16 three-phase power socket.

Returns the on-state of the CEE16 power socket

get_door_status(*door: int*) → hvl_ccb.dev.supercube2015.constants.DoorStatus

Get the status of a safety fence door. See `constants.DoorStatus` for possible returned door statuses.

Parameters `door` – the door number (1..3)

Returns the door status

get_earthing_manual(*number: int*) → bool

Get the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

Parameters `number` – number of the earthing stick (1..6)

Returns earthing stick manual status

get_earthing_status(*number: int*) → int

Get the status of an earthing stick, whether it is closed, open or undefined (moving).

Parameters `number` – number of the earthing stick (1..6)

Returns earthing stick status; see `constants.EarthingStickStatus`

get_measurement_ratio(*channel: int*) → float

Get the set measurement ratio of an AC/DC analog input channel. Every input channel has a divider ratio assigned during setup of the Supercube system. This ratio can be read out.

Attention: Supercube 2015 does not have a separate ratio for every analog input. Therefore there is only one ratio for `channel = 1`.

Parameters `channel` – number of the input channel (1..4)

Returns the ratio

get_measurement_voltage(*channel: int*) → float

Get the measured voltage of an analog input channel. The voltage read out here is already scaled by the configured divider ratio.

Attention: In contrast to the *new* Supercube, the old one returns here the input voltage read at the ADC. It is not scaled by a factor.

Parameters `channel` – number of the input channel (1..4)

Returns measured voltage

get_status() → int

Get the safety circuit status of the Supercube.

Returns the safety status of the supercube's state machine; see *constants.SafetyStatus*.

get_support_input(*port: int, contact: int*) → bool

Get the state of a support socket input.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital input read state

get_support_output(*port: int, contact: int*) → bool

Get the state of a support socket output.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)

Returns digital output read state

get_t13_socket(*port: int*) → bool

Read the state of a SEV T13 power socket.

Parameters **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*

Returns on-state of the power socket

horn(*state: bool*) → None

Turns acoustic horn on or off.

Parameters **state** – Turns horn on (True) or off (False)

operate(*state: bool*) → None

Set operate state. If the state is RedReady, this will turn on the high voltage and close the safety switches.

Parameters **state** – set operate state

quit_error() → None

Quits errors that are active on the Supercube.

read(*node_id: str*)

Local wrapper for the OPC UA communication protocol read method.

Parameters **node_id** – the id of the node to read.

Returns the value of the variable

ready(*state: bool*) → None

Set ready state. Ready means locket safety circuit, red lamps, but high voltage still off.

Parameters **state** – set ready state

set_cee16_socket(*state: bool*) → None

Switch the IEC CEE16 three-phase power socket on or off.

Parameters **state** – desired on-state of the power socket

Raises **ValueError** – if state is not of type bool

set_earthing_manual(*number: int, manual: bool*) → None

Set the manual status of an earthing stick. If an earthing stick is set to manual, it is closed even if the system is in states RedReady or RedOperate.

Parameters

- **number** – number of the earthing stick (1..6)
- **manual** – earthing stick manual status (True or False)

set_remote_control(*state: bool*) → None

Enable or disable remote control for the Supercube. This will effectively display a message on the touch-screen HMI.

Parameters **state** – desired remote control state

set_support_output(*port: int, contact: int, state: bool*) → None

Set the state of a support output socket.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **state** – is the desired state of the support output

set_support_output_impulse(*port: int, contact: int, duration: float = 0.2, pos_pulse: bool = True*) → None

Issue an impulse of a certain duration on a support output contact. The polarity of the pulse (On-wait-Off or Off-wait-On) is specified by the pos_pulse argument.

This function is blocking.

Parameters

- **port** – is the socket number (1..6)
- **contact** – is the contact on the socket (1..2)
- **duration** – is the length of the impulse in seconds
- **pos_pulse** – is True, if the pulse shall be HIGH, False if it shall be LOW

set_t13_socket(*port: int, state: bool*) → None

Set the state of a SEV T13 power socket.

Parameters

- **port** – is the socket number, one of *constants.T13_SOCKET_PORTS*
- **state** – is the desired on-state of the socket

start() → None

Starts the device. Sets the root node for all OPC read and write commands to the Siemens PLC object node which holds all our relevant objects and variables.

stop() → None

Stop the Supercube device. Deactivates the remote control and closes the communication protocol.

write(*node_id, value*) → None

Local wrapper for the OPC UA communication protocol write method.

Parameters

- **node_id** – the id of the node to read

- **value** – the value to write to the variable

class SupercubeConfiguration(namespace_index: int = 7)
 Bases: object

Configuration dataclass for the Supercube devices.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

namespace_index: int = 7

Namespace of the OPC variables, typically this is 3 (coming from Siemens)

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

class SupercubeOpcUaCommunication(config)

Bases: *hvl_ccb.comm.opc.OpcUaCommunication*

Communication protocol specification for Supercube devices.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class SupercubeOpcUaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], endpoint_name: str, port: int = 4845, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5)

Bases: *hvl_ccb.comm.opc.OpcUaCommunicationConfig*

Communication protocol configuration for OPC UA, specifications for the Supercube devices.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 4845

Port of the OPC UA server to connect to.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler =

<hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object>

Subscription handler for data change events

class SupercubeSubscriptionHandler

Bases: *hvl_ccb.comm.opc.OpcUaSubHandler*

OPC Subscription handler for datachange events and normal events specifically implemented for the Supercube devices.


datachange_notification(*node: opcua.common.node.Node*, *val*, *data*)

In addition to the standard operation (debug logging entry of the datachange), alarms are logged at INFO level using the alarm text.

Parameters

- **node** – the node object that triggered the datachange event
- **val** – the new value
- **data** –

hvl_ccb.dev.supercube2015.constants

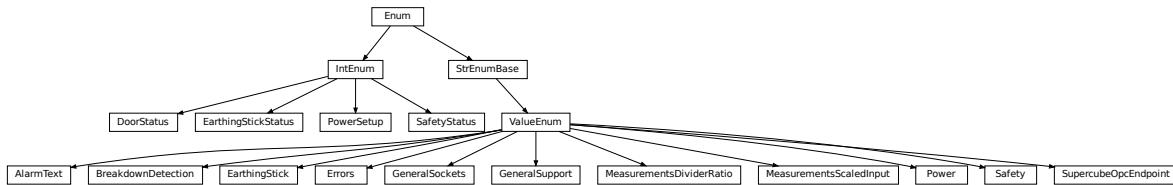

Constants, variable names for the Supercube OPC-connected devices.

```
class AlarmText(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum
```

This enumeration contains textual representations for all error classes (stop, warning and message) of the Supercube system. Use the `AlarmText.get()` method to retrieve the enum of an alarm number.

```
Alarm0 = 'No Alarm.'

Alarm1 = 'STOP Safety switch 1 error'
Alarm10 = 'STOP Earthing stick 2 error'
Alarm11 = 'STOP Earthing stick 3 error'
Alarm12 = 'STOP Earthing stick 4 error'
Alarm13 = 'STOP Earthing stick 5 error'
Alarm14 = 'STOP Earthing stick 6 error'
Alarm17 = 'STOP Source switch error'
Alarm19 = 'STOP Fence 1 error'
Alarm2 = 'STOP Safety switch 2 error'
Alarm20 = 'STOP Fence 2 error'
Alarm21 = 'STOP Control error'
Alarm22 = 'STOP Power outage'
Alarm3 = 'STOP Emergency Stop 1'
Alarm4 = 'STOP Emergency Stop 2'
Alarm5 = 'STOP Emergency Stop 3'
Alarm6 = 'STOP Door 1 lock supervision'
Alarm7 = 'STOP Door 2 lock supervision'
Alarm8 = 'STOP Door 3 lock supervision'
Alarm9 = 'STOP Earthing stick 1 error'

classmethod get(alarm: int)
```

Get the attribute of this enum for an alarm number.

Parameters `alarm` – the alarm number

Returns the enum for the desired alarm number

```
not_defined = 'NO ALARM TEXT DEFINED'
```

```
class BreakdownDetection(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Node ID strings for the breakdown detection.

activated = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.Breakdowndetection.connect'
Boolean read-only variable indicating whether breakdown detection and fast switchoff is enabled in the
system or not.

reset = 'hvl-ipc.WINAC.Support60utA'
Boolean writable variable to reset the fast switch-off. Toggle to re-enable.

triggered = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.Breakdowndetection.triggered'
Boolean read-only variable telling whether the fast switch-off has triggered. This can also be seen using
the safety circuit state, therefore no method is implemented to read this out directly.

class DoorStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum

Possible status values for doors.

closed = 2
Door is closed, but not locked.

error = 4
Door has an error or was opened in locked state (either with emergency stop or from the inside).

inactive = 0
not enabled in Supercube HMI setup, this door is not supervised.

locked = 3
Door is closed and locked (safe state).

open = 1
Door is open.

class EarthingStick(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.ValueEnum

Variable NodeID strings for all earthing stick statuses (read-only integer) and writable booleans for setting the
earthing in manual mode.

classmethod manual(number: int)
Get the manual enum attribute for an earthing stick number.

Parameters number – the earthing stick (1..6)

Returns the manual enum

manual_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.MANUAL'
manual_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.MANUAL'
manual_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.MANUAL'
manual_4 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.MANUAL'
manual_5 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.MANUAL'
manual_6 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.MANUAL'
status_1_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.CLOSE'
status_1_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.CONNECT'
```

```

status_1_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_1.OPEN'
status_2_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.CLOSE'
status_2_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.CONNECT'
status_2_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_2.OPEN'
status_3_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.CLOSE'
status_3_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.CONNECT'
status_3_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_3.OPEN'
status_4_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.CLOSE'
status_4_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.CONNECT'
status_4_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_4.OPEN'
status_5_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.CLOSE'
status_5_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.CONNECT'
status_5_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_5.OPEN'
status_6_closed = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.CLOSE'
status_6_connected = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.CONNECT'
status_6_open = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.STICK_6.OPEN'

```

classmethod status_closed(number: int)

Get the status enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status enum

classmethod status_connected(number: int)

Get the status enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status enum

classmethod status_open(number: int)

Get the status enum attribute for an earthing stick number.

Parameters **number** – the earthing stick (1..6)

Returns the status enum

```
class EarthingStickStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
```

Bases: aenum.IntEnum

Status of an earthing stick. These are the possible values in the status integer e.g. in EarthingStick.status_1.

closed = 1

Earthing is closed (safe).

error = 3

Earthing is in error, e.g. when the stick did not close correctly or could not open.

inactive = 0

Earthing stick is deselected and not enabled in safety circuit. To get out of this state, the earthing has to be enabled in the Supercube HMI setup.

open = 2

Earthing is open (not safe).

class Errors(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

Variable NodeID strings for information regarding error, warning and message handling.

quit = 'hvl-ipc.WINAC.SYSTEMSTATE.Faultconfirmation'

Writable boolean for the error quit button.

stop = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'

Boolean read-only variable telling if a stop is active.

stop_number = 'hvl-ipc.WINAC.SYSTEMSTATE.Errornumber'

class GeneralSockets(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

NodeID strings for the power sockets (3x T13 and 1xCEE16).

cee16 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.CEE16'

CEE16 socket (writeable boolean).

t13_1 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_1'

SEV T13 socket No. 1 (writable boolean).

t13_2 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_2'

SEV T13 socket No. 2 (writable boolean).

t13_3 = 'hvl-ipc.WINAC.SYSTEM_COMPONENTS.T13_3'

SEV T13 socket No. 3 (writable boolean).

class GeneralSupport(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: *hvl_ccb.utils.enum.ValueEnum*

NodeID strings for the support inputs and outputs.

in_1_1 = 'hvl-ipc.WINAC.Support1InA'

in_1_2 = 'hvl-ipc.WINAC.Support1InB'

in_2_1 = 'hvl-ipc.WINAC.Support2InA'

in_2_2 = 'hvl-ipc.WINAC.Support2InB'

in_3_1 = 'hvl-ipc.WINAC.Support3InA'

in_3_2 = 'hvl-ipc.WINAC.Support3InB'

in_4_1 = 'hvl-ipc.WINAC.Support4InA'

in_4_2 = 'hvl-ipc.WINAC.Support4InB'

in_5_1 = 'hvl-ipc.WINAC.Support5InA'

in_5_2 = 'hvl-ipc.WINAC.Support5InB'

in_6_1 = 'hvl-ipc.WINAC.Support6InA'

in_6_2 = 'hvl-ipc.WINAC.Support6InB'

classmethod **input**(*port, contact*)

Get the NodeID string for a support input.

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string

```
out_1_1 = 'hvl-ipc.WINAC.Support1OutA'
out_1_2 = 'hvl-ipc.WINAC.Support1OutB'
out_2_1 = 'hvl-ipc.WINAC.Support2OutA'
out_2_2 = 'hvl-ipc.WINAC.Support2OutB'
out_3_1 = 'hvl-ipc.WINAC.Support3OutA'
out_3_2 = 'hvl-ipc.WINAC.Support3OutB'
out_4_1 = 'hvl-ipc.WINAC.Support4OutA'
out_4_2 = 'hvl-ipc.WINAC.Support4OutB'
out_5_1 = 'hvl-ipc.WINAC.Support5OutA'
out_5_2 = 'hvl-ipc.WINAC.Support5OutB'
out_6_1 = 'hvl-ipc.WINAC.Support6OutA'
out_6_2 = 'hvl-ipc.WINAC.Support6OutB'
```

classmethod **output**(*port, contact*)

Get the NodeID string for a support output.

Parameters

- **port** – the desired port (1..6)
- **contact** – the desired contact at the port (1..2)

Returns the node id string

```
class MeasurementsDividerRatio(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: *hvl_ccb.utils.enum.ValueEnum*

Variable NodeID strings for the measurement input scaling ratios. These ratios are defined in the Supercube HMI setup and are provided in the python module here to be able to read them out, allowing further calculations.

classmethod **get**(*channel: int*)

Get the attribute for an input number.

Parameters **channel** – the channel number (1..4)

Returns the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.DivididerRatio'
```

```
class MeasurementsScaledInput(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: *hvl_ccb.utils.enum.ValueEnum*

Variable NodeID strings for the four analog BNC inputs for measuring voltage. The voltage returned in these variables is already scaled with the set ratio, which can be read using the variables in [MeasurementsDividerRatio](#).

classmethod get(channel: int)

Get the attribute for an input number.

Parameters **channel** – the channel number (1..4)

Returns the enum for the desired channel.

```
input_1 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI1Volt'
input_2 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI2Volt'
input_3 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI3Volt'
input_4 = 'hvl-ipc.WINAC.SYSTEM_INTERN.AI4Volt'
```

class Power(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: [hvl_ccb.utils.enum.ValueEnum](#)

Variable NodeID strings concerning power data.

current_primary = 'hvl-ipc.WINAC.SYSTEM_INTERN.FUCurrentprim'

Primary current in ampere, measured by the frequency converter. (read-only)

frequency = 'hvl-ipc.WINAC.FU.Frequency'

Frequency converter output frequency. (read-only)

setup = 'hvl-ipc.WINAC.FU.TrafoSetup'

Power setup that is configured using the Supercube HMI. The value corresponds to the ones in [PowerSetup](#). (read-only)

voltage_max = 'hvl-ipc.WINAC.FU.maxVoltagekV'

Maximum voltage allowed by the current experimental setup. (read-only)

voltage_primary = 'hvl-ipc.WINAC.SYSTEM_INTERN.FUVoltageprim'

Primary voltage in volts, measured by the frequency converter at its output. (read-only)

voltage_slope = 'hvl-ipc.WINAC.FU.dUdt_-1'

Voltage slope in V/s.

voltage_target = 'hvl-ipc.WINAC.FU.SOLL'

Target voltage setpoint in V.

class PowerSetup(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: [aenum.IntEnum](#)

Possible power setups corresponding to the value of variable [Power.setup](#).

AC_DoubleStage_150kV = 3

AC voltage with two MWB transformers, one at 100kV and the other at 50kV, resulting in a total maximum voltage of 150kV.

AC_DoubleStage_200kV = 4

AC voltage with two MWB transformers both at 100kV, resulting in a total maximum voltage of 200kV

AC_SingleStage_100kV = 2

AC voltage with MWB transformer set to 100kV maximum voltage.

AC_SingleStage_50kV = 1

AC voltage with MWB transformer set to 50kV maximum voltage.

DC_DoubleStage_280kV = 7

DC voltage with two AC transformers set to 100kV AC each, resulting in 280kV DC in total (or a single stage transformer with Greinacher voltage doubling rectifier)

DC_SingleStage_140kV = 6

DC voltage with one AC transformer set to 100kV AC, resulting in 140kV DC

External = 0

External power supply fed through blue CEE32 input using isolation transformer and safety switches of the Supercube, or using an external safety switch attached to the Supercube Type B.

Internal = 5

Internal usage of the frequency converter, controlling to the primary voltage output of the supercube itself (no measurement transformer used)

class Safety(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: [hvl_ccb.utils.enum.ValueEnum](#)

NodeID strings for the basic safety circuit status and green/red switches “ready” and “operate”.

horn = 'hvl-ipc.WINAC.SYSTEM_INTERN.hornen'

Writeable boolean to manually turn on or off the horn

status_error = 'hvl-ipc.WINAC.SYSTEMSTATE.ERROR'**status_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN'****status_ready_for_red = 'hvl-ipc.WINAC.SYSTEMSTATE.ReadyForRed'**

Status is a read-only integer containing the state number of the supercube-internal state machine. The values correspond to numbers in [SafetyStatus](#).

status_red = 'hvl-ipc.WINAC.SYSTEMSTATE.RED'**switchto_green = 'hvl-ipc.WINAC.SYSTEMSTATE.GREEN_REQUEST'****switchto_operate = 'hvl-ipc.WINAC.SYSTEMSTATE.switchon'**

Writable boolean for switching to Red Operate (locket, HV on) state.

switchto_ready = 'hvl-ipc.WINAC.SYSTEMSTATE.RED_REQUEST'

Writable boolean for switching to Red Ready (locked, HV off) state.

class SafetyStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: [aenum.IntEnum](#)

Safety status values that are possible states returned from [hvl_ccb.dev.supercube.base.Supercube.get_status\(\)](#). These values correspond to the states of the Supercube’s safety circuit statemachine.

Error = 6

System is in error mode.

GreenNotReady = 1

System is safe, lamps are green and some safety elements are not in place such that it cannot be switched to red currently.

GreenReady = 2

System is safe and all safety elements are in place to be able to switch to *ready*.

Initializing = 0

System is initializing or booting.

QuickStop = 5

Fast turn off triggered and switched off the system. Reset FSO to go back to a normal state.

RedOperate = 4

System is locked in red state and in *operate* mode, i.e. high voltage on.

RedReady = 3

System is locked in red state and *ready* to go to *operate* mode.

```
class SupercubeOpcEndpoint(value=<no_arg>, names=None, module=None, type=None, start=1,
                             boundary=None)
```

Bases: *hvl_ccb.utils.enum.ValueEnum*

OPC Server Endpoint strings for the supercube variants.


A = 'OPC.SimaticNET.S7'

B = 'OPC.SimaticNET.S7'

T13_SOCKET_PORTS = (1, 2, 3)

Port numbers of SEV T13 power socket

[hvl_ccb.dev.supercube2015.typ_a](#)

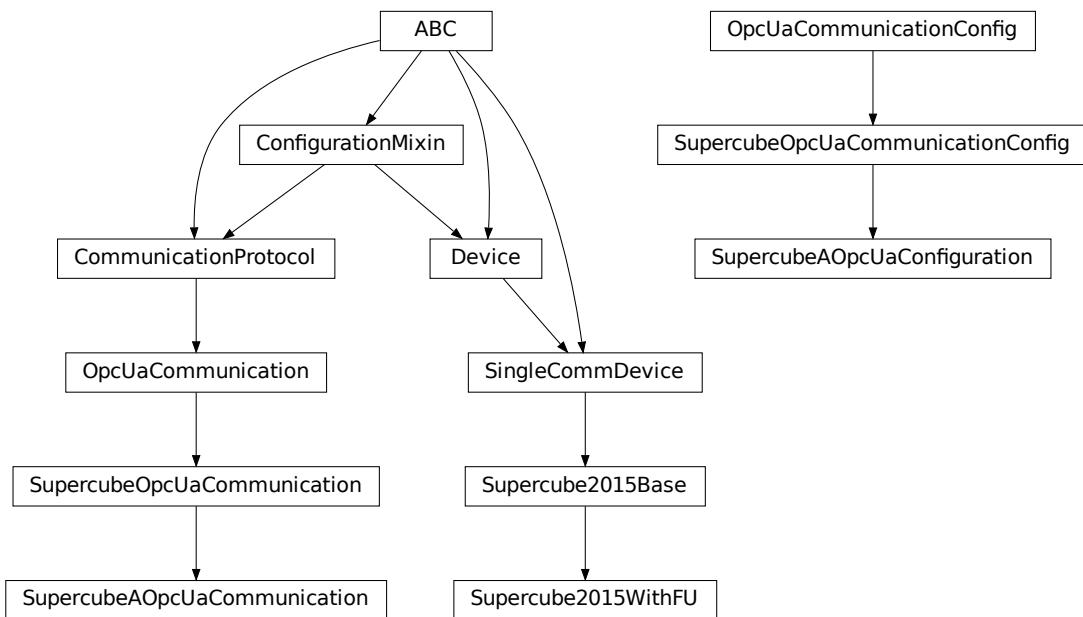

Supercube Typ A module.

```
class Supercube2015WithFU(com, dev_config=None)
```

Bases: *hvl_ccb.dev.supercube2015.base.Supercube2015Base*

Variant A of the Supercube with frequency converter.

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

fso_reset() → None
 Reset the fast switch off circuitry to go back into normal state and allow to re-enable operate mode.

get_frequency() → float
 Read the electrical frequency of the current Supercube setup.
Returns the frequency in Hz

get_fso_active() → bool
 Get the state of the fast switch off functionality. Returns True if it is enabled, False otherwise.
Returns state of the FSO functionality

get_max_voltage() → float
 Reads the maximum voltage of the setup and returns in V.
Returns the maximum voltage of the setup in V.

get_power_setup() → *hvl_ccb.dev.supercube2015.constants.PowerSetup*
 Return the power setup selected in the Supercube's settings.
Returns the power setup

get_primary_current() → float
 Read the current primary current at the output of the frequency converter (before transformer).
Returns primary current in A

get_primary_voltage() → float
 Read the current primary voltage at the output of the frequency converter (before transformer).
Returns primary voltage in V

get_target_voltage() → float
 Gets the current setpoint of the output voltage value in V. This is not a measured value but is the corresponding function to *set_target_voltage()*.
Returns the setpoint voltage in V.

set_slope(*slope: float*) → None
 Sets the dV/dt slope of the Supercube frequency converter to a new value in V/s.
Parameters **slope** – voltage slope in V/s (0..15'000)

set_target_voltage(*volt_v: float*) → None
 Set the output voltage to a defined value in V.
Parameters **volt_v** – the desired voltage in V

class SupercubeA0pcUaCommunication(*config*)
 Bases: *hvl_ccb.dev.supercube2015.base.SupercubeOpcUaCommunication*

static config_cls()
 Return the default configdataclass class.
Returns a reference to the default configdataclass class

class SupercubeA0pcUaConfiguration(*host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], endpoint_name: str = 'OPC.SimaticNET.S7', port: int = 4845, sub_handler: hvl_ccb.comm.opc.OpcUaSubHandler = <hvl_ccb.dev.supercube2015.base.SupercubeSubscriptionHandler object at 0x7fb049bc3dd0>, update_period: int = 500, wait_timeout_retry_sec: Union[int, float] = 1, max_timeout_retry_nr: int = 5*)
 Bases: *hvl_ccb.dev.supercube2015.base.SupercubeOpcUaCommunicationConfig*

```
endpoint_name: str = 'OPC.SimaticNET.S7'  
Endpoint of the OPC server, this is a path like 'OPCUA/SimulationServer'  
force_value(fieldname, value)  
Forces a value to a dataclass field despite the class being frozen.  
NOTE: you can define post_force_value method with same signature as this method to do extra processing  
after value has been forced on fieldname.  
Parameters

- fieldname – name of the field
- value – value to assign

classmethod keys() → Sequence[str]  
Returns a list of all configdataclass fields key-names.  
Returns a list of strings containing all keys.  
classmethod optional_defaults() → Dict[str, object]  
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified  
on instantiation.  
Returns a list of strings containing all optional keys.  
classmethod required_keys() → Sequence[str]  
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on  
instantiation.  
Returns a list of strings containing all required keys.
```


Module contents

Supercube package with implementation for the old system version from 2015 based on Siemens WinAC soft-PLC on
an industrial 32bit Windows computer.

`hvl_ccb.dev.technix`

Submodules

`hvl_ccb.dev.technix.base`

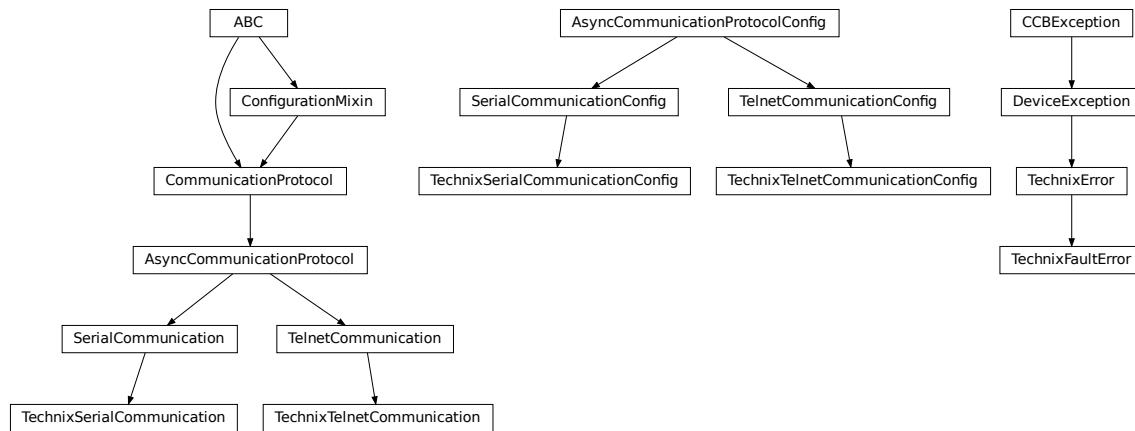

Communication and auxiliary classes for Technix

`exception TechnixError`

Bases: `hvl_ccb.dev.base.DeviceException`

Technix related errors.

`exception TechnixFaultError`

Bases: `hvl_ccb.dev.technix.base.TechinxError`

Raised when the fault flag was detected while the interlock is closed

`class TechnixSerialCommunication(configuration)`

Bases: `hvl_ccb.dev.technix.base._TechnixCommunication, hvl_ccb.comm.serial.SerialCommunication`

Serial communication for Technix

`static config_cls()`

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class TechnixSerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, port:
                                         Optional[str] = None, baudrate: int = 9600, parity: Union[str,
                                         hvl_ccb.comm.serial.SerialCommunicationParity] =
                                         SerialCommunicationParity.NONE, stopbits: Union[int, float,
                                         hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                         SerialCommunicationStopbits.ONE, bytesize: Union[int,
                                         hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                         SerialCommunicationBytesize.EIGHTBITS, timeout: Union[int,
                                         float] = 2)
```

Bases: `hvl_ccb.dev.technix.base._TechnixCommunicationConfig`, `hvl_ccb.comm.serial.SerialCommunicationConfig`

Configuration for the serial communication for Technix

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

class TechnixTelnetCommunication(*configuration*)

Bases: `hvl_ccb.comm.telnet.TelnetCommunication`, `hvl_ccb.dev.technix.base._TechnixCommunication`

Telnet communication for Technix

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

```
class TechnixTelnetCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',
                                         encoding_error_handling: str = 'strict',
                                         wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                         default_n_attempts_read_text_nonempty: int = 10, host:
                                         Optional[Union[str, ipaddress.IPv4Address,
                                         ipaddress.IPv6Address]] = None, port: int = 4660, timeout:
                                         Union[int, float] = 0.2)
Bases:      hvl_ccb.dev.technix.base._TechnixCommunicationConfig,  hvl_ccb.comm.telnet.
              TelnetCommunicationConfig
```

Configuration for the telnet communication for Technix

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

port: int = 4660


Port at which Technix is listening

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

`hvl_ccb.dev.technix.device`

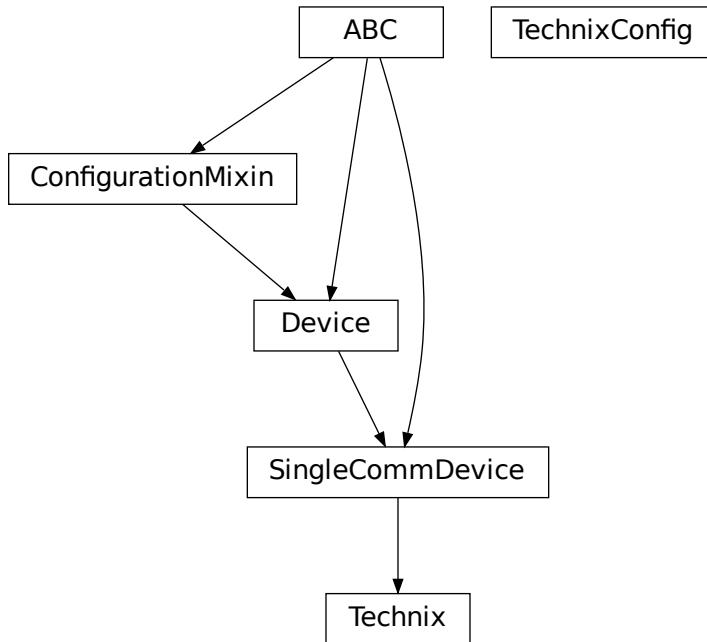

its corresponding configuration class

```

class Technix(com, dev_config)
  Bases: hvl_ccb.dev.base.SingleCommDevice
  Device class to control capacitor chargers from Technix

  static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

  property current: Optional[Union[int, float]]
    Actual current of the output in A

  default_com_cls() → Union[Type[hvl_ccb.dev.technix.base.TechnixSerialCommunication],
    Type[hvl_ccb.dev.technix.base.TechnixTelnetCommunication]]
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

  property inhibit: Optional[bool]
    Is the output of the voltage inhibited? The output stage can still be active.

  property is_started: bool
    Is the device started?

  property max_current: Union[int, float]
    Maximal output current of the hardware in A
  
```

The device class *Technix* and

```

property max_voltage: Union[int, float]
    Maximal output voltage of the hardware in V

property open_interlock: Optional[bool]
    Is the interlock open? (in safe mode)

property output: Optional[bool]
    State of the high voltage output

query_status(*, _retry: bool = False)
    Query the status of the device.

Returns This function returns nothing

property remote: Optional[bool]
    Is the device in remote control mode?

start()
    Start the device and set it into the remote controllable mode. The high voltage is turn off, and the status poller is started.

property status: Optional[hvl_ccb.dev.technix.base._Status]
    The status of the device with the different states as sub-fields

stop()
    Stop the device. The status poller is stopped and the high voltage output is turn off.

property voltage: Optional[Union[int, float]]
    Actual voltage at the output in V

property voltage_regulation: Optional[bool]
    Status if the output is in voltage regulation mode (or current regulation)

class TechnixConfig(communication_channel:
    Union[Type[hvl_ccb.dev.technix.base.TechnixSerialCommunication],
    Type[hvl_ccb.dev.technix.base.TechnixTelnetCommunication]], max_voltage: Union[int,
    float], max_current: Union[int, float], polling_interval_sec: Union[int, float] = 4,
    post_stop_pause_sec: Union[int, float] = 1, register_pulse_time: Union[int, float] = 0.1,
    read_output_while_polling: bool = False)

Bases: object

clean_values()
    Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

communication_channel:
Union[Type[hvl_ccb.dev.technix.base.TechnixSerialCommunication],
Type[hvl_ccb.dev.technix.base.TechnixTelnetCommunication]]
    communication channel between computer and Technix

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing after value has been forced on fieldname.

Parameters

- fieldname – name of the field
- value – value to assign

is_configdataclass = True

```

```
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

max_current: Union[int, float]
    Maximal Output current

max_voltage: Union[int, float]
    Maximal Output voltage

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

polling_interval_sec: Union[int, float] = 4
    Polling interval in s to maintain to watchdog of the device

post_stop_pause_sec: Union[int, float] = 1
    Time to wait after stopping the device

read_output_while_polling: bool = False
    Read output voltage and current within the polling event

register_pulse_time: Union[int, float] = 0.1
    Time for pulsing a register

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.
```


Module contents

Device classes for “RS 232” and “Ethernet” interfaces, which are used to control power supplies from Technix. Manufacturer homepage: <https://www.technix-hv.com>

The regulated power supplies series and capacitor chargers series from Technix are series of low and high voltage direct current power supplies as well as capacitor chargers. The class *Technix* is tested with a CCR10KV-7,5KJ via an ethernet connection as well as a CCR15-P-2500-OP via a serial connection. Check the code carefully before using it with other devices or device series

This Python package may support the following interfaces from Technix:

- Remote Interface RS232
- Ethernet Remote Interface
- Optic Fiber Remote Interface

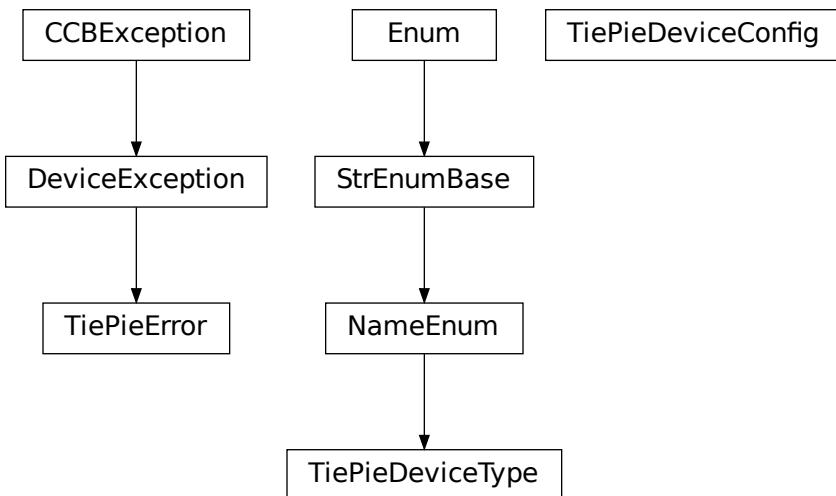hvl_ccb.dev.tiepie**Submodules****hvl_ccb.dev.tiepie.base**

```
class TiePieDeviceConfig(serial_number: int, require_block_measurement_support: bool = True,
                        n_max_try_get_device: int = 10, wait_sec_retry_get_device: Union[int, float] =
                        1.0, is_data_ready_polling_interval_sec: Union[int, float] = 0.01)
```

Bases: `object`

Configuration dataclass for TiePie

`clean_values()`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- `fieldname` – name of the field
- `value` – value to assign

`is_configdataclass = True`

`is_data_ready_polling_interval_sec: Union[int, float] = 0.01`

`classmethod keys() → Sequence[str]`

Returns a list of all configdataclass fields key-names.

`Returns` a list of strings containing all keys.

```
n_max_try_get_device: int = 10
classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

require_block_measurement_support: bool = True
classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

serial_number: int
wait_sec_retry_get_device: Union[int, float] = 1.0
class TiePieDeviceType(value=<no_arg>, names=None, module=None, type=None, start=1,
                       boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
TiePie device type.

GENERATOR = 2
I2C = 4
OSCILLOSCOPE = 1
exception TiePieError
Bases: hvl_ccb.dev.base.DeviceException
Error of the class TiePie

get_device_by_serial_number(serial_number: int, device_type: Union[str, Tuple[int,
                           hvl_ccb.dev.tiepie.base._LtpDeviceReturnType]], n_max_try_get_device: int =
                           10, wait_sec_retry_get_device: float = 1.0) →
                           hvl_ccb.dev.tiepie.base._LtpDeviceReturnType
Open and return handle of TiePie device with a given serial number
```

Parameters

- **serial_number** – int serial number of the device
- **device_type** – a *TiePieDeviceType* instance containing device identifier (int number) and its corresponding class, both from *libtiepie*, or a string name of such instance
- **n_max_try_get_device** – maximal number of device list updates (int number)
- **wait_sec_retry_get_device** – waiting time in seconds between retries (int number)

Returns Instance of a *libtiepie* device class according to the specified *device_type*

Raises

- **TiePieError** – when there is no device with given serial number
- **ValueError** – when *device_type* is not an instance of *TiePieDeviceType*

wrap_libtiepie_exception(func: Callable) → Callable


Decorator wrapper for *libtiepie* methods that use *libtiepie.library.check_last_status_raise_on_error()* calls.

Parameters **func** – Function or method to be wrapped

Raises `TiePieError` – instead of `LibTiePieException` or one of its subtypes.

Returns whatever `func` returns

hvl_ccb.dev.tiepie.channel

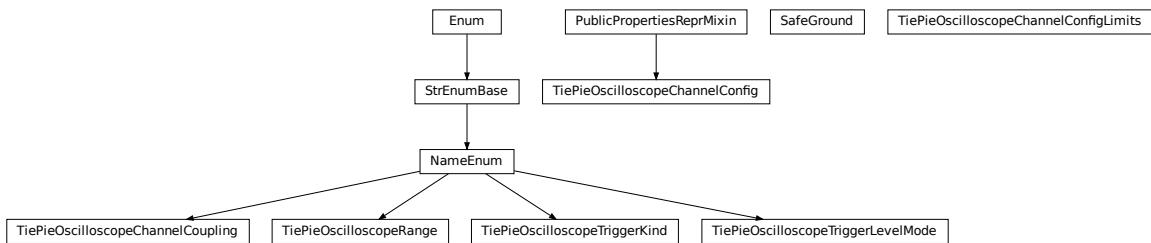

class SafeGround

Bases: `object`

Class that dynamically adds the `safe_ground_enabled` attribute getter/setter if the bound oscilloscope has the safe ground option.

```

class TiePieOscilloscopeChannelConfig(ch_number: int, channel:
    libtiepie.oscilloscopechannel.OscilloscopeChannel)
Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin

Oscilloscope's channel configuration, with cleaning of values in properties setters as well as setting and reading them on and from the device's channel.

static clean_coupling(coupling: Union[str,
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling]) →
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling

static clean_enabled(enabled: bool) → bool

clean_input_range(input_range: Union[float, hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeRange]) →
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeRange

clean_probe_offset(probe_offset: float) → float

static clean_trigger_enabled(trigger_enabled)

clean_trigger_hysteresis(trigger_hysteresis: float) → float

static clean_trigger_kind(trigger_kind: Union[str,
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind]) →
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind

clean_trigger_level(trigger_level: Union[int, float]) → float

static clean_trigger_level_mode(level_mode: Union[str,
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode]) →
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode

property coupling: hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelCoupling
property enabled: bool
  
```

```
property has_safe_ground: bool
    Check whether bound oscilloscope device has “safe ground” option

    Returns bool: 1=safe ground available

property input_range: hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeRange
property probe_offset: float
property trigger_enabled: bool
property trigger_hysteresis: float
property trigger_kind: hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerKind
property trigger_level: float
property trigger_level_mode:
    hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeTriggerLevelMode

class TiePieOscilloscopeChannelConfigLimits(osc_channel:
                                             libtiepie.oscilloscopechannel.OscilloscopeChannel)
Bases: object

Default limits for oscilloscope channel parameters.

class TiePieOscilloscopeChannelCoupling(value=<no_arg>, names=None, module=None, type=None,
                                         start=1, boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ACA = 8
ACV = 2
DCA = 4
DCV = 1

class TiePieOscilloscopeRange(value=<no_arg>, names=None, module=None, type=None, start=1,
                               boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

EIGHTY_VOLT = 80
EIGHT_HUNDRED_MILLI_VOLT = 0.8
EIGHT_VOLT = 8
FORTY_VOLT = 40
FOUR_HUNDRED_MILLI_VOLT = 0.4
FOUR_VOLT = 4
TWENTY_VOLT = 20
TWO_HUNDRED_MILLI_VOLT = 0.2
TWO_VOLT = 2

static suitable_range(value)
```

```
class TiePieOscilloscopeTriggerKind(value=<no_arg>, names=None, module=None, type=None, start=1,
                                     boundary=None)
```

Bases: `hvl_ccb.utils.enum.NameEnum`

An enumeration.

ANY = 16

FALLING = 2

RISING = 1

RISING_OR_FALLING = 16

```
class TiePieOscilloscopeTriggerLevelMode(value=<no_arg>, names=None, module=None, type=None,
                                         start=1, boundary=None)
```

Bases: `hvl_ccb.utils.enum.NameEnum`


An enumeration.

ABSOLUTE = 2

RELATIVE = 1

UNKNOWN = 0

[hvl_ccb.dev.tiepie.device](#)

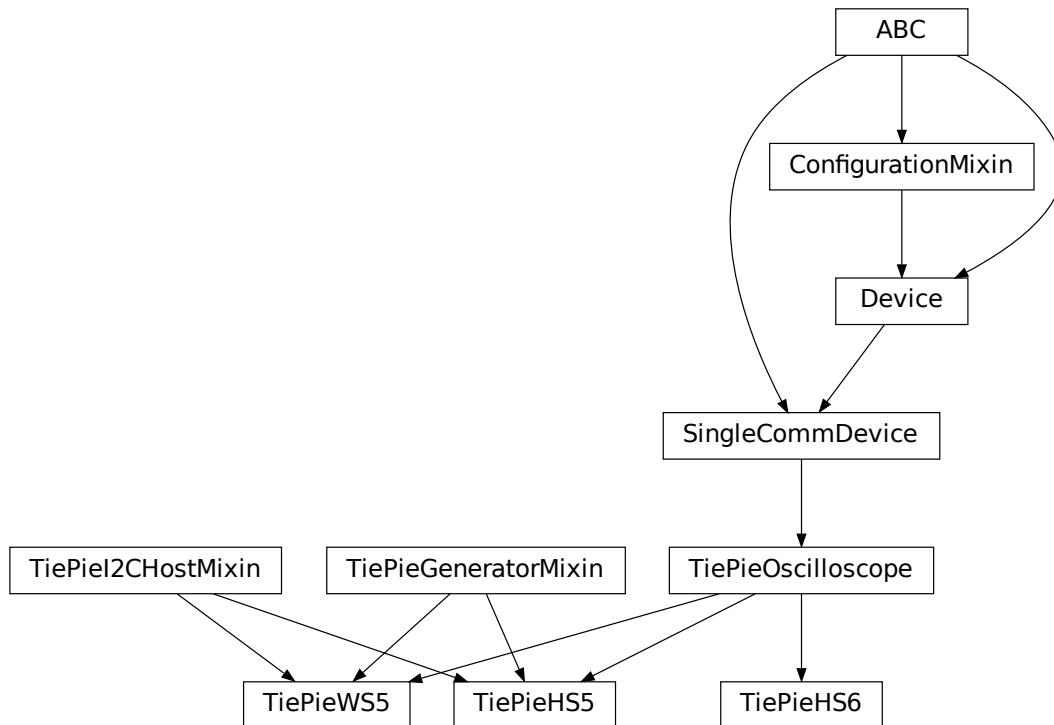

TiePie devices.

```
class TiePieHS5(com, dev_config)
Bases:      hvl_ccb.dev.tiepie.i2c.TiePieI2CHostMixin,    hvl_ccb.dev.tiepie.generator.
TiePieGeneratorMixin, hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope

TiePie HS5 device.


class TiePieHS6(com, dev_config)
Bases: hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope

TiePie HS6 DIFF device.

class TiePieWS5(com, dev_config)
Bases:      hvl_ccb.dev.tiepie.i2c.TiePieI2CHostMixin,    hvl_ccb.dev.tiepie.generator.
TiePieGeneratorMixin, hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscope

TiePie WS5 device.
```

hvl_ccb.dev.tiepie.generator

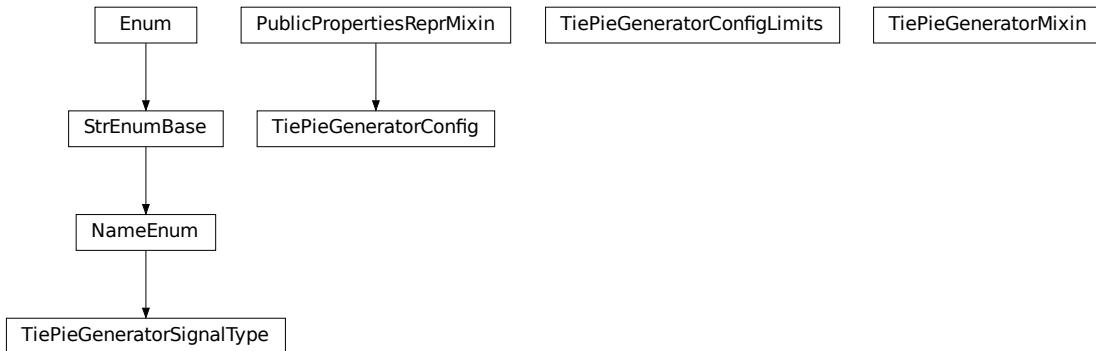

```
class TiePieGeneratorConfig(dev_gen: libtiepie.generator.Generator)
Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin

Generator's configuration with cleaning of values in properties setters.

property amplitude: float
clean_amplitude(amplitude: float) → float
static clean_enabled(enabled: bool) → bool
clean_frequency(frequency: float) → float
clean_offset(offset: float) → float
static clean_signal_type(signal_type: Union[int,
                                              hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType]) →
                                              hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType
clean_waveform(waveform: numpy.ndarray[Any, numpy.dtype[numpy.typing._generic_alias.ScalarType]]) →
                                              numpy.ndarray[Any, numpy.dtype[numpy.typing._generic_alias.ScalarType]]
property enabled: bool
```

```

property frequency: float
property offset: float
property signal_type: hvl_ccb.dev.tiepie.generator.TiePieGeneratorSignalType
property waveform: Optional[numpy.ndarray[Any,
numpy.dtype[numpy.typing._generic_alias.ScalarType]]]

class TiePieGeneratorConfigLimits(dev_gen: libtiepie.generator.Generator)
Bases: object

Default limits for generator parameters.

class TiePieGeneratorMixin(com, dev_config)
Bases: object

TiePie Generator sub-device.

A wrapper for the libtiepie.generator.Generator class. To be mixed in with TiePieOscilloscope base class.

config_gen: Optional[hvl_ccb.dev.tiepie.generator.TiePieGeneratorConfig]
Generator's dynamical configuration.

generator_start()
Start signal generation.

generator_stop()
Stop signal generation.

start() → None
Start the Generator.

stop() → None
Stop the generator.


class TiePieGeneratorSignalType(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ARBITRARY = 32
DC = 8
NOISE = 16
PULSE = 64
SINE = 1
SQUARE = 4
TRIANGLE = 2
UNKNOWN = 0

```

`hvl_ccb.dev.tiepie.i2c`

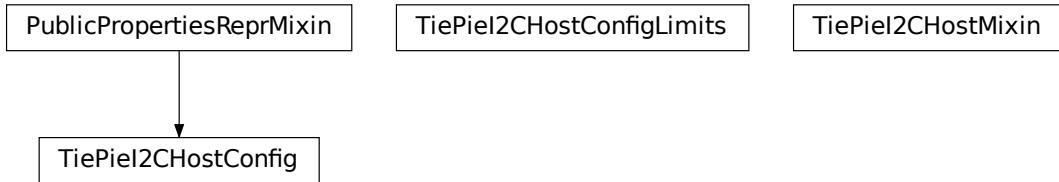

```
class TiePieI2CHostConfig(dev_i2c: libtiepie.i2chost.I2CHost)
    Bases: hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin
    I2C Host's configuration with cleaning of values in properties setters.

class TiePieI2CHostConfigLimits(dev_i2c: libtiepie.i2chost.I2CHost)
    Bases: object
    Default limits for I2C host parameters.


class TiePieI2CHostMixin(com, dev_config)
    Bases: object
    TiePie I2CHost sub-device.
    A wrapper for the libtiepie.i2chost.I2CHost class. To be mixed in with TiePieOscilloscope base class.

config_i2c: Optional[hvl_ccb.dev.tiepie.i2c.TiePieI2CHostConfig]
    I2C host's dynamical configuration.

start() → None
    Start the I2C Host.

stop() → None
    Stop the I2C host.
```

hvl_ccb.dev.tiepie.oscilloscope

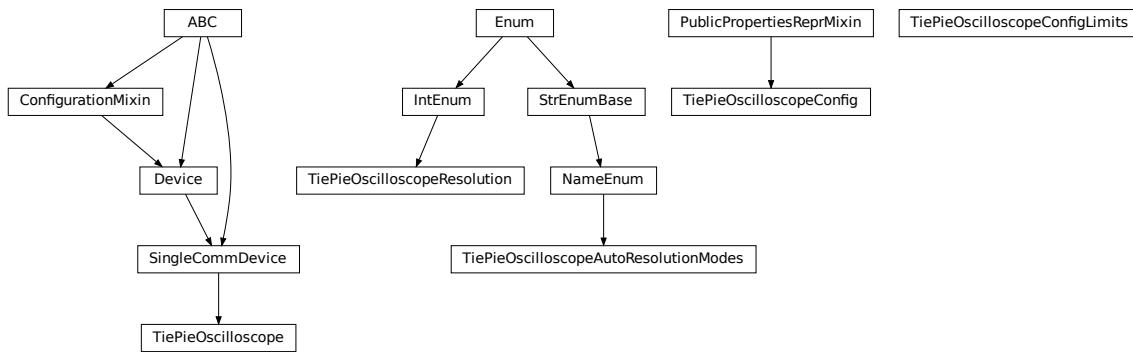

```
class TiePieOscilloscope(com, dev_config)
Bases: hvl\_ccb.dev.base.SingleCommDevice
```

TiePie oscilloscope.

A wrapper for TiePie oscilloscopes, based on the class *libtiepie.oscilloscope.Oscilloscope* with simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and its channels, including extra validation and typing hints support for configurations.

Note that, in contrast to *libtiepie* library, since all physical TiePie devices include an oscilloscope, this is the base class for all physical TiePie devices. The additional TiePie sub-devices: “Generator” and “I2CHost”, are mixed-in to this base class in subclasses.

The channels use *I..N* numbering (not *0..N-1*), as in, e.g., the Multi Channel software.

property channels_enabled: Generator[int, None, None]

Yield numbers of enabled channels.

Returns Numbers of enabled channels

collect_measurement_data(timeout: Optional[Union[int, float]] = 0) → Optional[numpy.ndarray[Any, numpy.dtype[numpy.typing._generic_alias.ScalarType]]]

Try to collect the data from TiePie; return *None* if data is not ready.

Parameters **timeout** – The timeout to wait until data is available. This option makes this function blocking the code. *timeout = None* blocks the code infinitely till data will be available. Per default, the *timeout* is set to *0*: The function will not block.

Returns Measurement data of only enabled channels and time vector in a 2D-*numpy.ndarray* with float sample data; or *None* if there is no data available.

static config_cls() → Type[[hvl_ccb.dev.tiepie.base.TiePieDeviceConfig](#)]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

config_osc: Optional[[hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeConfig](#)]

Oscilloscope’s dynamical configuration.

config_osc_channel_dict: Dict[int,

[hvl_ccb.dev.tiepie.channel.TiePieOscilloscopeChannelConfig](#)

Channel configuration. A *dict* mapping actual channel number, numbered *I..N*, to channel configuration.

The channel info is dynamically read from the device only on the first `start()`; beforehand the `dict` is empty.

static default_com_cls() → Type[*hvl_ccb.comm.base.NullCommunicationProtocol*]

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

force_trigger() → None

Forces the TiePie to trigger with a software sided trigger event.

Return None

Raises *TiePieError* – when device is not started or status of underlying device gives an error

is_measurement_data_ready() → bool

Reports if TiePie has data which is ready to collect

Returns if the data is ready to collect.

Raises *TiePieError* – when device is not started or status of underlying device gives an error

is_triggered() → bool

Reports if TiePie has triggered. Maybe data is not yet available. One can check with the function `is_measurement_data_ready()`.

Returns if a trigger event occurred

static list_devices() → libtiepie.devicelist.DeviceList

List available TiePie devices.

Returns libtiepie up to date list of devices

property n_channels

Number of channels in the oscilloscope.

Returns Number of channels.

start() → None

Start the oscilloscope.

start_measurement() → None

Start a measurement using set configuration.

Raises *TiePieError* – when device is not started or status of underlying device gives an error

stop() → None

Stop the oscilloscope.

class TiePieOscilloscopeAutoResolutionModes(*value=<no_arg>*, *names=None*, *module=None*,
type=None, *start=1*, *boundary=None*)

Bases: *hvl_ccb.utils.enum.NameEnum*

An enumeration.

ALL = 4

DISABLED = 1

NATIVEONLY = 2

UNKNOWN = 0

class TiePieOscilloscopeConfig(*dev_osc: libtiepie.oscilloscope.Oscilloscope*)

Bases: *hvl_ccb.dev.tiepie.utils.PublicPropertiesReprMixin*

Oscilloscope's configuration with cleaning of values in properties setters.

```

property auto_resolution_mode:
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes

static clean_auto_resolution_mode(auto_resolution_mode: Union[int,
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes]) →
    →
        hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeAutoResolutionModes

clean_pre_sample_ratio(pre_sample_ratio: float) → float
clean_record_length(record_length: Union[int, float]) → int
static clean_resolution(resolution: Union[int,
    hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution]) →
        hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution

clean_sample_frequency(sample_frequency: float) → float
clean_trigger_timeout(trigger_timeout: Optional[Union[int, float]]) → float
property pre_sample_ratio: float
property record_length: int
property resolution: hvl_ccb.dev.tiepie.oscilloscope.TiePieOscilloscopeResolution
property sample_frequency: float
property trigger_timeout: Optional[float]

class TiePieOscilloscopeConfigLimits(dev_osc: libtiepie.oscilloscope.Oscilloscope)
Bases: object

Default limits for oscilloscope parameters.

class TiePieOscilloscopeResolution(value=<no_arg>, names=None, module=None, type=None, start=1,
    boundary=None)
Bases: aenum.IntEnum

An enumeration.

EIGHT_BIT = 8
FOURTEEN_BIT = 14
SIXTEEN_BIT = 16
TWELVE_BIT = 12

```

[hvl_ccb.dev.tiepie.utils](#)

PublicPropertiesReprMixin

```

class PublicPropertiesReprMixin
Bases: object

```

General purpose utility mixin that overwrites object representation to a one analogous to *dataclass* instances, but using public properties and their values instead of *fields*.

Module contents

This module is a wrapper around LibTiePie SDK devices; see <https://www.tiepie.com/en/libtiepie-sdk> .

The device classes adds simplifications for starting of the device (using serial number) and managing mutable configuration of both the device and oscilloscope's channels. This includes extra validation and typing hints support.

Extra installation

LibTiePie SDK library is available only on Windows and on Linux.

To use this LibTiePie SDK devices wrapper:

1. install the hvl_ccb package with a tiepie extra feature:

```
$ pip install "hvl_ccb[tiepie]"
```

this will install the Python bindings for the library.

2. install the library

- on Linux: follow instructions in <https://www.tiepie.com/en/libtiepie-sdk/linux> ;
- on Windows: the additional DLL is included in Python bindings package.

Troubleshooting

On a Windows system, if you encounter an OSError like this:

```
...
self._handle = _dlopen(self._name, mode)
OSError: [WinError 126] The specified module could not be found
```


most likely the python-libtiepie package was installed in your site-packages/ directory as a python-libtiepie-*.egg file via python setup.py install or python setup.py develop command. In such case uninstall the library and re-install it using pip:

```
$ pip uninstall python-libtiepie
$ pip install python-libtiepie
```

This should create libtiepie/ folder. Alternatively, manually move the folder libtiepie/ from inside of the .egg archive file to the containing it site-packages/ directory (PyCharm's Project tool window supports reading and extracting from .egg archives).

Submodules

`hvl_ccb.dev.base`

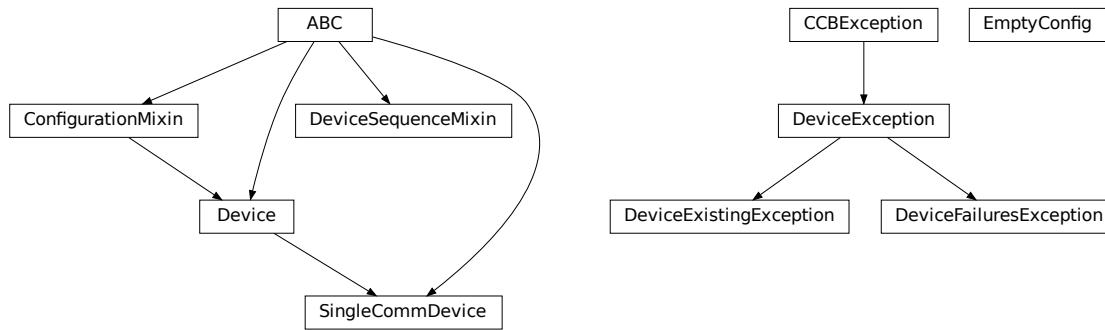

Module with base classes for devices.

class Device(`dev_config=None`)
Bases: `hvl_ccb.configuration.ConfigurationMixin, abc.ABC`

Base class for devices. Implement this class for a concrete device, such as measurement equipment or voltage sources.

Specifies the methods to implement for a device.

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

abstract start() → None
Start or restart this Device. To be implemented in the subclass.

abstract stop() → None
Stop this Device. To be implemented in the subclass.

exception DeviceException
Bases: `hvl_ccb.exception.CCBException`

exception DeviceExistingException
Bases: `hvl_ccb.dev.base.DeviceException`

Exception to indicate that a device with that name already exists.

exception DeviceFailuresException(`failures: Dict[str, Exception], *args`)
Bases: `hvl_ccb.dev.base.DeviceException`

Exception to indicate that one or several devices failed.

failures: Dict[str, Exception]
A dictionary of named devices failures (exceptions).

class DeviceSequenceMixin(`devices: Dict[str, hvl_ccb.dev.base.Device]`)
Bases: `abc.ABC`

Mixin that can be used on a device or other classes to provide facilities for handling multiple devices in a sequence.

add_device(*name: str, device: hvl_ccb.dev.base.Device*) → None

Add a new device to the device sequence.

Parameters

- **name** – is the name of the device.
- **device** – is the instantiated Device object.

Raises *DeviceExistingException* –

devices_failed_start: Dict[str, hvl_ccb.dev.base.Device]

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices_failed_stop*.

devices_failed_stop: Dict[str, hvl_ccb.dev.base.Device]

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices_failed_start*.

get_device(*name: str*) → hvl_ccb.dev.base.Device

Get a device by name.

Parameters **name** – is the name of the device.

Returns the device object from this sequence.

get_devices() → List[Tuple[str, hvl_ccb.dev.base.Device]]

Get list of name, device pairs according to current sequence.

Returns A list of tuples with name and device each.

remove_device(*name: str*) → hvl_ccb.dev.base.Device

Remove a device from this sequence and return the device object.

Parameters **name** – is the name of the device.

Returns device object or *None* if such device was not in the sequence.

Raises **ValueError** – when device with given name was not found

start() → None

Start all devices in this sequence in their added order.

Raises *DeviceFailuresException* – if one or several devices failed to start

stop() → None

Stop all devices in this sequence in their reverse order.

Raises *DeviceFailuresException* – if one or several devices failed to stop

class EmptyConfig

Bases: object

Empty configuration dataclass that is the default configuration for a Device.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

class SingleCommDevice(com, dev_config=None)

Bases: *hvl_ccb.dev.base.Device*, abc.ABC

Base class for devices with a single communication protocol.

property com

Get the communication protocol of this device.

Returns an instance of CommunicationProtocol subtype

abstract static default_com_cls() → Type[*hvl_ccb.comm.base.CommunicationProtocol*]

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device


start() → None

Open the associated communication protocol.

stop() → None

Close the associated communication protocol.

hvl_ccb.dev.crylas

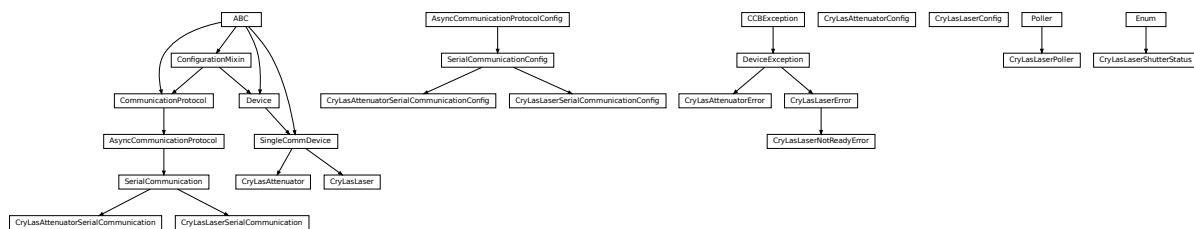

Device classes for a CryLas pulsed laser controller and a CryLas laser attenuator, using serial communication.

There are three modes of operation for the laser 1. Laser-internal hardware trigger (default): fixed to 20 Hz and max energy per pulse. 2. Laser-internal software trigger (for diagnosis only). 3. External trigger: required for arbitrary pulse energy or repetition rate. Switch to “external” on the front panel of laser controller for using option 3.

After switching on the laser with `laser_on()`, the system must stabilize for some minutes. Do not apply abrupt changes of pulse energy or repetition rate.

Manufacturer homepage: https://www.crylas.de/products/pulsed_laser.html

class CryLasAttenuator(*com, dev_config=None*)
Bases: `hvl_ccb.dev.base.SingleCommDevice`

Device class for the CryLas laser attenuator.

property attenuation: `Union[int, float]`

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

set_attenuation(*percent: Union[int, float]*) → None

Set the percentage of attenuated light (inverse of set_transmission). :param percent: percentage of attenuation, number between 0 and 100 :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

set_init_attenuation()

Sets the attenuation to its configured initial/default value

Raises `SerialCommunicationIOError` – when communication port is not opened

set_transmission(*percent: Union[int, float]*) → None

Set the percentage of transmitted light (inverse of set_attenuation). :param percent: percentage of transmitted light :raises ValueError: if param percent not between 0 and 100 :raises SerialCommunicationIOError: when communication port is not opened :raises CryLasAttenuatorError: if the device does not confirm success

start() → None

Open the com, apply the config value ‘init_attenuation’

Raises `SerialCommunicationIOError` – when communication port cannot be opened

property transmission: `Union[int, float]`

class CryLasAttenuatorConfig(*init_attenuation: Union[int, float] = 0, response_sleep_time: Union[int, float] = 1*)
Bases: `object`

Device configuration dataclass for CryLas attenuator.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

```

init_attenuation: Union[int, float] = 0
is_configdataclass = True

classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

response_sleep_time: Union[int, float] = 1

exception CryLasAttenuatorError
    Bases: hvl_ccb.dev.base.DeviceException
    General error with the CryLas Attenuator.

class CryLasAttenuatorSerialCommunication(configuration)
    Bases: hvl_ccb.comm.serial.SerialCommunication
    Specific communication protocol implementation for the CryLas attenuator. Already predefines device-specific protocol parameters in config.

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

class CryLasAttenuatorSerialCommunicationConfig(terminator: bytes = b'', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)
    Bases: hvl_ccb.comm.serial.SerialCommunicationConfig

baudrate: int = 9600
    Baudrate for CryLas attenuator is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
    One byte is eight bits long

```

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, *hvl_ccb.comm.serial.SerialCommunicationParity*] = 'N'

CryLas attenuator does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, *hvl_ccb.comm.serial.SerialCommunicationStopbits*] = 1

CryLas attenuator uses one stop bit

terminator: bytes = b''

No terminator

timeout: Union[int, float] = 3

use 3 seconds timeout as default

class CryLasLaser(*com*, *dev_config*=None)

Bases: *hvl_ccb.dev.base.SingleCommDevice*

CryLas laser controller device class.

class AnswersShutter(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *aenum.Enum*

Standard answers of the CryLas laser controller to 'Shutter' command passed via *com*.

CLOSED = 'Shutter inaktiv'

OPENED = 'Shutter aktiv'

class AnswersStatus(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: *aenum.Enum*

Standard answers of the CryLas laser controller to 'STATUS' command passed via *com*.

ACTIVE = 'STATUS: Laser active'

HEAD = 'STATUS: Head ok'

```

INACTIVE = 'STATUS: Laser inactive'
READY = 'STATUS: System ready'
TEC1 = 'STATUS: TEC1 Regulation ok'
TEC2 = 'STATUS: TEC2 Regulation ok'

class LaserStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: aenum.Enum

    Status of the CryLas laser

    READY_ACTIVE = 2
    READY_INACTIVE = 1
    UNREADY_INACTIVE = 0
    property is_inactive
    property is_ready

class RepetitionRates(value=<no_arg>, names=None, module=None, type=None, start=1,
                      boundary=None)
    Bases: aenum.IntEnum

    Repetition rates for the internal software trigger in Hz

    HARDWARE = 0
    SOFTWARE_INTERNAL_SIXTY = 60
    SOFTWARE_INTERNAL_TEN = 10
    SOFTWARE_INTERNAL_TWENTY = 20

ShutterStatus
    alias of hvl\_ccb.dev.crylas.CryLasLaserShutterStatus

close_shutter() → None
    Close the laser shutter.

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • CryLasLaserError – if success is not confirmed by the device

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

get_pulse_energy_and_rate() → Tuple[int, int]
    Use the debug mode, return the measured pulse energy and rate.

    Returns (energy in micro joule, rate in Hz)

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • CryLasLaserError – if the device does not answer the query

```

laser_off() → None

Turn the laser off.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

laser_on() → None

Turn the laser on.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserNotReadyError* – if the laser is not ready to be turned on
- *CryLasLaserError* – if success is not confirmed by the device

open_shutter() → None

Open the laser shutter.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

set_init_shutter_status() → None

Open or close the shutter, to match the configured shutter_status.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

set_pulse_energy(energy: int) → None

Sets the energy of pulses (works only with external hardware trigger). Proceed with small energy steps, or the regulation may fail.

Parameters `energy` – energy in micro joule

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if the device does not confirm success

set_repetition_rate(rate: Union[int, hvl_ccb.dev.crylas.CryLasLaser.RepetitionRates]) → None

Sets the repetition rate of the internal software trigger.

Parameters `rate` – frequency (Hz) as an integer

Raises

- *ValueError* – if rate is not an accepted value in RepetitionRates Enum
- *SerialCommunicationIOError* – when communication port is not opened
- *CryLasLaserError* – if success is not confirmed by the device

start() → None

Opens the communication protocol and configures the device.

Raises *SerialCommunicationIOError* – when communication port cannot be opened

stop() → None

Stops the device and closes the communication protocol.

Raises

- **SerialCommunicationIOError** – if com port is closed unexpectedly
- **CryLasLaserError** – if laser_off() or close_shutter() fail

property target_pulse_energy**update_laser_status()** → NoneUpdate the laser status to *LaserStatus.NOT_READY* or *LaserStatus.INACTIVE* or *LaserStatus.ACTIVE*.

Note: laser never explicitly says that it is not ready (*LaserStatus.NOT_READY*) in response to ‘STATUS’ command. It only says that it is ready (heated-up and implicitly inactive/off) or active (on). If it’s not either of these then the answer is *Answers.HEAD*. Moreover, the only time the laser explicitly says that its status is inactive (*Answers.INACTIVE*) is after issuing a ‘LASER OFF’ command.

Raises **SerialCommunicationIOError** – when communication port is not opened

update_repetition_rate() → None

Query the laser repetition rate.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

update_shutter_status() → None

Update the shutter status (OPENED or CLOSED)

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

update_target_pulse_energy() → None

Query the laser pulse energy.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **CryLasLaserError** – if success is not confirmed by the device

wait_until_ready() → None

Block execution until the laser is ready

Raises **CryLasLaserError** – if the polling thread stops before the laser is ready

```
class CryLasLaserConfig(calibration_factor: Union[int, float] = 4.35, polling_period: Union[int, float] = 12,
                        polling_timeout: Union[int, float] = 300, auto_laser_on: bool = True,
                        init_shutter_status: Union[int, hvl_ccb.dev.crylas.CryLasLaserShutterStatus] =
                        CryLasLaserShutterStatus.CLOSED)
```

Bases: **object**

Device configuration dataclass for the CryLas laser controller.

ShutterStatusalias of *hvl_ccb.dev.crylas.CryLasLaserShutterStatus***auto_laser_on:** bool = True**calibration_factor:** Union[int, float] = 4.35

`clean_values()`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

`init_shutter_status: Union[int, hvl_ccb.dev.crylas.CryLasLaserShutterStatus] = 0`

`is_configdataclass = True`

`classmethod keys() → Sequence[str]`

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

`classmethod optional_defaults() → Dict[str, object]`

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

`polling_period: Union[int, float] = 12`

`polling_timeout: Union[int, float] = 300`

`classmethod required_keys() → Sequence[str]`

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception CryLasLaserError

Bases: `hvl_ccb.dev.base.DeviceException`

General error with the CryLas Laser.

exception CryLasLaserNotReadyError

Bases: `hvl_ccb.dev.crylas.CryLasLaserError`

Error when trying to turn on the CryLas Laser before it is ready.

class CryLasLaserPoller(spoll_handler: Callable, check_handler: Callable, check_laser_status_handler:

`Callable, polling_delay_sec: Union[int, float] = 0, polling_interval_sec: Union[int, float] = 1, polling_timeout_sec: Optional[Union[int, float]] = None)`

Bases: `hvl_ccb.dev.utils.Poller`

Poller class for polling the laser status until the laser is ready.

Raises

- `CryLasLaserError` – if the timeout is reached before the laser is ready
- `SerialCommunicationIOError` – when communication port is closed.

class CryLasLaserSerialCommunication(configuration)

Bases: `hvl_ccb.comm.serial.SerialCommunication`

Specific communication protocol implementation for the CryLas laser controller. Already predefines device-specific protocol parameters in config.

READ_TEXT_SKIP_PREFIXES = ('>', 'MODE:')

Prefixes of lines that are skipped when read from the serial port.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(cmd: str, prefix: str, post_cmd: Optional[str] = None) → str

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns the line in question.

Parameters

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer
- **post_cmd** – optional additional command to send after the query

Returns line in question as a string

Raises *SerialCommunicationIOError* – when communication port is not opened

query_all(cmd: str, prefix: str)

Send a command, then read the com until a line starting with prefix, or an empty line, is found. Returns a list of successive lines starting with prefix.

Parameters

- **cmd** – query message to send to the device
- **prefix** – start of the line to look for in the device answer

Returns line in question as a string

Raises *SerialCommunicationIOError* – when communication port is not opened

read() → str

Read first line of text from the serial port that does not start with any of *self.READ_TEXT_SKIP_PREFIXES*.

Returns String read from the serial port; '' if there was nothing to read.

Raises *SerialCommunicationIOError* – when communication port is not opened

class CryLasLaserSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 19200, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 10)

Bases: *hvl_ccb.comm.serial.SerialCommunicationConfig*

```
baudrate: int = 19200
    Baudrate for CryLas laser is 19200 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
    One byte is eight bits long

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing
    after value has been forced on fieldname.
```

Parameters

- **fieldname** – name of the field
- **value** – value to assign

```
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.
```

```
classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified
    on instantiation.

    Returns a list of strings containing all optional keys.
```

```
parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'
    CryLas laser does not use parity
```

```
classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on
    instantiation.

    Returns a list of strings containing all required keys.
```

```
stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
    CryLas laser uses one stop bit
```


```
terminator: bytes = b'\n'
    The terminator is LF
```

```
timeout: Union[int, float] = 10
    use 10 seconds timeout as default (a long timeout is needed!)
```

```
class CryLasLaserShutterStatus(value=<no_arg>, names=None, module=None, type=None, start=1,
                                 boundary=None)
Bases: aenum.Enum

Status of the CryLas laser shutter

CLOSED = 0
OPENED = 1
```

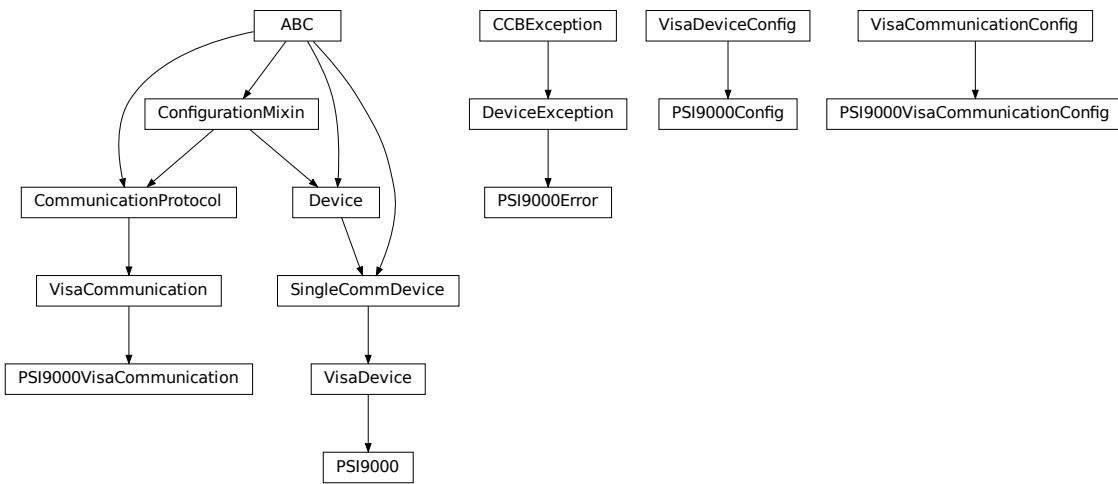hvl_ccb.dev.ea_psi9000

Device class for controlling a Elektro Automatik PSI 9000 power supply over VISA.

It is necessary that a backend for pyvisa is installed. This can be NI-Visa oder pyvisa-py (up to now, all the testing was done with NI-Visa)

```

class PSI9000(com: Union[hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunication,
hvl_ccb.dev.ea_psi9000.PSI9000VisaCommunicationConfig, dict], dev_config:
Optional[Union[hvl_ccb.dev.ea_psi9000.PSI9000Config, dict]] = None)
  
```

Bases: `hvl_ccb.dev.visa.VisaDevice`

Elektro Automatik PSI 9000 power supply.

MS_NOMINAL_CURRENT = 2040

MS_NOMINAL_VOLTAGE = 80

SHUTDOWN_CURRENT_LIMIT = 0.1

SHUTDOWN_VOLTAGE_LIMIT = 0.1

check_master_slave_config() → None

Checks if the master / slave configuration and initializes if successful

Returns `PSI9000Error` – if master-slave configuration failed

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Return the default communication protocol for this device type, which is `VisaCommunication`.

Returns the `VisaCommunication` class

get_output() → bool

Reads the current state of the DC output of the source. Returns True, if it is enabled, false otherwise.

Returns the state of the DC output

get_system_lock() → bool

Get the current lock state of the system. The lock state is true, if the remote control is active and false, if not.

Returns the current lock state of the device

get_ui_lower_limits() → Tuple[float, float]

Get the lower voltage and current limits. A lower power limit does not exist.

Returns Umin in V, Imin in A

get_uip_upper_limits() → Tuple[float, float, float]

Get the upper voltage, current and power limits.

Returns Umax in V, Imax in A, Pmax in W

get_voltage_current_setpoint() → Tuple[float, float]

Get the voltage and current setpoint of the current source.

Returns Uset in V, Iset in A

measure_voltage_current() → Tuple[float, float]

Measure the DC output voltage and current

Returns Umeas in V, Imeas in A

set_lower_limits(voltage_limit: Optional[float] = None, current_limit: Optional[float] = None) → None

Set the lower limits for voltage and current. After writing the values a check is performed if the values are set correctly.

Parameters

- **voltage_limit** – is the lower voltage limit in V
- **current_limit** – is the lower current limit in A

Raises `PSI9000Error` – if the limits are out of range

set_output(target_onstate: bool) → None

Enables / disables the DC output.

Parameters **target_onstate** – enable or disable the output power

Raises `PSI9000Error` – if operation was not successful

set_system_lock(lock: bool) → None

Lock / unlock the device, after locking the control is limited to this class unlocking only possible when voltage and current are below the defined limits

Parameters **lock** – True: locking, False: unlocking

set_upper_limits(voltage_limit: Optional[float] = None, current_limit: Optional[float] = None, power_limit: Optional[float] = None) → None

Set the upper limits for voltage, current and power. After writing the values a check is performed if the values are set. If a parameter is left blank, the maximum configurable limit is set.

Parameters

- **voltage_limit** – is the voltage limit in V
- **current_limit** – is the current limit in A
- **power_limit** – is the power limit in W

Raises `PSI9000Error` – if limits are out of range

set_voltage_current(*volt*: float, *current*: float) → None

Set voltage and current setpoints.

After setting voltage and current, a check is performed if writing was successful.

Parameters

- **volt** – is the setpoint voltage: 0.81.6 V (1.02 * 0-80 V) (absolute max, can be smaller if limits are set)
- **current** – is the setpoint current: 0..2080.8 A (1.02 * 0 - 2040 A) (absolute max, can be smaller if limits are set)

Raises `PSI9000Error` – if the desired setpoint is out of limits

start() → None

Start this device.

stop() → None

Stop this device. Turns off output and lock, if enabled.

class PSI9000Config(*spoll_interval*: Union[int, float] = 0.5, *spoll_start_delay*: Union[int, float] = 2, *power_limit*: Union[int, float] = 43500, *voltage_lower_limit*: Union[int, float] = 0.0, *voltage_upper_limit*: Union[int, float] = 10.0, *current_lower_limit*: Union[int, float] = 0.0, *current_upper_limit*: Union[int, float] = 2040.0, *wait_sec_system_lock*: Union[int, float] = 0.5, *wait_sec_settings_effect*: Union[int, float] = 1, *wait_sec_initialisation*: Union[int, float] = 2)

Bases: `hvl_ccb.dev.visa.VisaDeviceConfig`

Elektro Automatik PSI 9000 power supply device class. The device is communicating over a VISA TCP socket.

Using this power supply, DC voltage and current can be supplied to a load with up to 2040 A and 80 V (using all four available units in parallel). The maximum power is limited by the grid, being at 43.5 kW available through the CEE63 power socket.

clean_values() → None

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

current_lower_limit: Union[int, float] = 0.0

Lower current limit in A, depending on the experimental setup.

current_upper_limit: Union[int, float] = 2040.0

Upper current limit in A, depending on the experimental setup.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

power_limit: Union[int, float] = 43500

Power limit in W depending on the experimental setup. With 3x63A, this is 43.5kW. Do not change this value, if you do not know what you are doing. There is no lower power limit.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

voltage_lower_limit: Union[int, float] = 0.0

Lower voltage limit in V, depending on the experimental setup.

voltage_upper_limit: Union[int, float] = 10.0

Upper voltage limit in V, depending on the experimental setup.

wait_sec_initialisation: Union[int, float] = 2

wait_sec_settings_effect: Union[int, float] = 1

wait_sec_system_lock: Union[int, float] = 0.5

exception PSI9000Error

Bases: *hvl_ccb.dev.base.DeviceException*

Base error class regarding problems with the PSI 9000 supply.

class PSI9000VisaCommunication(configuration)

Bases: *hvl_ccb.comm.visa.VisaCommunication*

Communication protocol used with the PSI 9000 power supply.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class PSI9000VisaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], interface_type: Union[str, hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType] = InterfaceType.TCPIP_SOCKET, board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = '')

Bases: *hvl_ccb.comm.visa.VisaCommunicationConfig*

Visa communication protocol config dataclass with specification for the PSI 9000 power supply.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

interface_type: Union[str, *hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType*]
= 1

Interface type of the VISA connection, being one of InterfaceType.

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.


Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

hvl_ccb.dev.fug

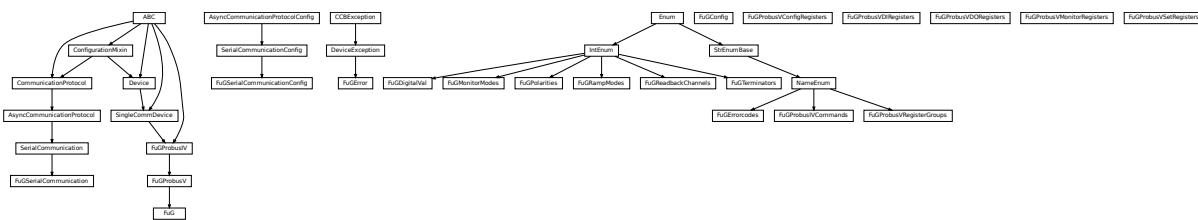

Device classes for “Probus V - ADDAT30” Interfaces which are used to control power supplies from FuG Elektronik GmbH

This interface is used for many FuG power units. Manufacturer homepage: <https://www.fug-elektronik.de>

The Professional Series of Power Supplies from FuG is a series of low, medium and high voltage direct current power supplies as well as capacitor chargers. The class FuG is tested with a HCK 800-20 000 in Standard Mode. The addressable mode is not implemented. Check the code carefully before using it with other devices. Manufacturer homepage: <https://www.fug-elektronik.de/netzgeraete/professional-series/>

The documentation of the interface from the manufacturer can be found here: https://www.fug-elektronik.de/wp-content/uploads/download/de/SOFTWARE/Probus_V.zip

The provided classes support the basic and some advanced commands. The commands for calibrating the power supplies are not implemented, as they are only for very special purposes and should not be used by “normal” customers.

class FuG(com, dev_config=None)

Bases: *hvl_ccb.dev.fug.FuGProbusV*

FuG power supply device class.

The power supply is controlled over a FuG ADDA Interface with the PROBUS V protocol

property config_status: *hvl_ccb.dev.fug.FuGProbusVConfigRegisters*

Returns the registers for the registers with the configuration and status values

Returns FuGProbusVConfigRegisters

property current: *hvl_ccb.dev.fug.FuGProbusVSetRegisters*

Returns the registers for the current output

Returns

property current_monitor: *hvl_ccb.dev.fug.FuGProbusVMonitorRegisters*

Returns the registers for the current monitor.

A typically usage will be “self.current_monitor.value” to measure the output current

Returns

property di: *hvl_ccb.dev.fug.FuGProbusVDIRegisters*

Returns the registers for the digital inputs

Returns FuGProbusVDIRegisters

identify_device() → None

Identify the device nominal voltage and current based on its model number.

Raises *SerialCommunicationIOError* – when communication port is not opened

property max_current: *Union[int, float]*

Returns the maximal current which could provided within the test setup

Returns

property max_current_hardware: *Union[int, float]*

Returns the maximal current which could provided with the power supply

Returns

property max_voltage: *Union[int, float]*

Returns the maximal voltage which could provided within the test setup

Returns

property max_voltage_hardware: *Union[int, float]*

Returns the maximal voltage which could provided with the power supply

Returns

property on: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the output switch to turn the output on or off

Returns FuGProbusVDORegisters

property outX0: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X0

Returns FuGProbusVDORegisters

property outX1: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X1

Returns FuGProbusVDORegisters

property outX2: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital output X2

Returns FuGProbusVDORegisters

property outXCMD: *hvl_ccb.dev.fug.FuGProbusVDORegisters*

Returns the registers for the digital outputX-CMD

Returns FuGProbusVDORegisters

start(*max_voltage=0, max_current=0*) → None
 Opens the communication protocol and configures the device.

Parameters

- **max_voltage** – Configure here the maximal permissible voltage which is allowed in the given experimental setup
- **max_current** – Configure here the maximal permissible current which is allowed in the given experimental setup

property voltage: *hvl_ccb.dev.fug.FuGProbusVSetRegisters*

Returns the registers for the voltage output

Returns

property voltage_monitor: *hvl_ccb.dev.fug.FuGProbusVMonitorRegisters*

Returns the registers for the voltage monitor.

A typically usage will be “self.voltage_monitor.value” to measure the output voltage

Returns

class FuGConfig(*wait_sec_stop_commands: Union[int, float] = 0.5*)

Bases: *object*

Device configuration dataclass for FuG power supplies.

clean_values()

force_value(*fieldname, value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

wait_sec_stop_commands: Union[int, float] = 0.5

Time to wait after subsequent commands during stop (in seconds)

```
class FuGDigitalVal(value)
    Bases: enum.IntEnum
    An enumeration.

    NO = 0
    OFF = 0
    ON = 1
    YES = 1

exception FuGError(*args, **kwargs)
    Bases: hvl_ccb.dev.base.DeviceException
    Error with the FuG voltage source.

    errorcode: str
        Errorcode from the Probus, see documentation of Probus V chapter 5. Errors with three-digit errorcodes
        are thrown by this python module.

class FuGErrorcodes(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl_ccb.utils.enum.NameEnum

    The power supply can return an errorcode. These errorcodes are handled by this class. The original errorcodes
    from the source are with one or two digits, see documentation of Probus V chapter 5. All three-digit errorcodes
    are from this python module.

    E0 = ('no error', 'standard response on each command')
    E1 = ('no data available', 'Customer tried to read from GPIB but there were no data
          prepared. (IBIG50 sent command ~T2 to ADDA)')
    E10 = ('unknown SCPI command', 'This SCPI command is not implemented')
    E100 = ('Command is not implemented', 'You tried to execute a command, which is not
            implemented or does not exist')
    E106 = ('The rampstate is a read-only register', 'You tried to write data to the
             register, which can only give you the status of the ramping.')
    E11 = ('not allowed Trigger-on-Talk', 'Not allowed attempt to Trigger-on-Talk (~T1)
           while ADDA was in addressable mode.')
    E115 = ('The given index to select a digital value is out of range', 'Only integer
             values between 0 and 1 are allowed.')
    E12 = ('invalid argument in ~Tn command', 'Only ~T1 and ~T2 is implemented.')
    E125 = ('The given index to select a ramp mode is out of range', 'Only integer
             values between 0 and 4 are allowed.')
    E13 = ('invalid N-value', 'Register > K8 contained an invalid value. Error code is
           output on an attempt to query data with ? or ~T1')
    E135 = ('The given index to select the readback channel is out of range', 'Only
             integer values between 0 and 6 are allowed.')
    E14 = ('register is write only', 'Some registers can only be writte to (i.e.> H0)')
    E145 = ('The given value for the AD-conversion is unknown', 'Valid values for the
             ad-conversion are integer values from "0" to "7".')
    E15 = ('string too long', 'i.e.serial number string too long during calibration')
```

```

E155 = ('The given value to select a polarity is out range.', 'The value should be 0 or 1.')
E16 = ('wrong checksum', 'checksum over command string was not correct, refer also to 4.4 of the Probus V documentation')
E165 = ('The given index to select the terminator string is out of range', '')
E2 = ('unknown register type', "No valid register type after '>'")
E206 = ('This status register is read-only', 'You tried to write data to this register, which can only give you the actual status of the corresponding digital output.')
E306 = ('The monitor register is read-only', 'You tried to write data to a monitor, which can only give you measured data.')
E4 = ('invalid argument', 'The argument of the command was rejected .i.e. malformed number')
E5 = ('argument out of range', 'i.e. setvalue higher than type value')
E504 = ('Empty string as response', 'The connection is broken.')
E505 = ('The returned register is not the requested.', 'Maybe the connection is overburden.')
E6 = ('register is read only', 'Some registers can only be read but not written to. (i.e. monitor registers)')
E666 = ('You cannot overwrite the most recent error in the interface of the power supply. But, well: You created an error anyway...', '')
E7 = ('Receive Overflow', 'Command string was longer than 50 characters.')
E8 = ('EEPROM is write protected', 'Write attempt to calibration data while the write protection switch was set to write protected.')
E9 = ('address error', 'A non addressed command was sent to ADDA while it was in addressable mode (and vice versa.)')
raise_()

class FuGMonitorModes(value)
Bases: enum.IntEnum
An enumeration.

T1MS = 1
    15 bit + sign, 1 ms integration time

T200MS = 6
    typ. 19 bit + sign, 200 ms integration time

T20MS = 3
    17 bit + sign, 20 ms integration time

T256US = 0
    14 bit + sign, 256 us integration time

T40MS = 4
    17 bit + sign, 40 ms integration time

T4MS = 2
    15 bit + sign, 4 ms integration time

```

```
T800MS = 7
typ. 20 bit + sign, 800 ms integration time

T80MS = 5
typ. 18 bit + sign, 80 ms integration time

class FuGPolarities(value)
Bases: enum.IntEnum

An enumeration.

NEGATIVE = 1
POSITIVE = 0

class FuGProbusIV(com, dev_config=None)
Bases: hvl_ccb.dev.base.SingleCommDevice, abc.ABC

FuG Probus IV device class

Sends basic SCPI commands and reads the answer. Only the special commands and PROBUS IV instruction set
is implemented.

command(command: hvl_ccb.dev.fug.FuGProbusIVCommands, value=None) → str

Parameters

- command – one of the commands given within FuGProbusIVCommands
- value – an optional value, depending on the command

Returns a String if a query was performed

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

output_off() → None
Switch DC voltage output off.

reset() → None
Reset of the interface: All setvalues are set to zero

abstract start()
Open the associated communication protocol.

stop() → None
Close the associated communication protocol.

class FuGProbusIVCommands(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum

An enumeration.

ADMODE = ('S', (<enum 'FuGMonitorModes'>, <class 'int'>))
CURRENT = ('I', (<class 'int'>, <class 'float'>))
EXECUTE = ('X', None)
```

```

EXECUTEONX = ('G', (<enum 'FuGDigitalVal'>, <class 'int'>))
    Wait for "X" to execute pending commands

ID = ('*IDN?', None)

OUTPUT = ('F', (<enum 'FuGDigitalVal'>, <class 'int'>))

POLARITY = ('P', (<enum 'FuGPolarities'>, <class 'int'>))

QUERY = ('?', None)

READBACKCHANNEL = ('N', (<enum 'FuGReadbackChannels'>, <class 'int'>))

RESET = ('=', None)

TERMINATOR = ('Y', (<enum 'FuGTerminators'>, <class 'int'>))

VOLTAGE = ('U', (<class 'int'>, <class 'float'>))

XOUTPUTS = ('R', <class 'int'>)
    TODO: the possible values are limited to 0..13

```

class FuGProbusV(*com, dev_config=None*)
Bases: *hvl_ccb.dev.fug.FuGProbusIV*

FuG Probus V class which uses register based commands to control the power supplies

get_register(*register: str*) → str
get the value from a register

Parameters **register** – the register from which the value is requested

Returns the value of the register as a String

set_register(*register: str, value: Union[int, float, str]*) → None
generic method to set value to register

Parameters

- **register** – the name of the register to set the value
- **value** – which should be written to the register

class FuGProbusVConfigRegisters(*fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups*)
Bases: object

Configuration and Status values, acc. 4.2.5

property execute_on_x: *hvl_ccb.dev.fug.FuGDigitalVal*
status of Execute-on-X

Returns FuGDigitalVal of the status

property most_recent_error: *hvl_ccb.dev.fug.FuGErrorcodes*
Reads the Error-Code of the most recent command

Return FuGError

Raises *FuGError* – if code is not “E0”

property readback_data: *hvl_ccb.dev.fug.FuGReadbackChannels*
Preselection of readout data for Trigger-on-Talk

Returns index for the readback channel

property srq_mask: *int*
SRQ-Mask, Service-Request Enable status bits for SRQ 0: no SRQ Bit 2: SRQ on change of status to CC
Bit 1: SRQ on change to CV

Returns representative integer value

property srq_status: str
SRQ-Statusbyte output as a decimal number: Bit 2: PS is in CC mode Bit 1: PS is in CV mode

Returns representative string

property status: str
Statusbyte as a string of 0/1. Combined status (compatibel to Probus IV), MSB first: Bit 7: I-REG Bit 6: V-REG Bit 5: ON-Status Bit 4: 3-Reg Bit 3: X-Stat (polarity) Bit 2: Cal-Mode Bit 1: unused Bit 0: SEL-D

Returns string of 0/1

property terminator: hvl_ccb.dev.fug.FuGTerminators
Terminator character for answer strings from ADDA

Returns FuGTerminators

class FuGProbusVDIRegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Digital Inputs acc. 4.2.4

property analog_control: hvl_ccb.dev.fug.FuGDigitalVal
Returns shows 1 if power supply is controlled by the analog interface

property calibration_mode: hvl_ccb.dev.fug.FuGDigitalVal
Returns shows 1 if power supply is in calibration mode

property cc_mode: hvl_ccb.dev.fug.FuGDigitalVal
Returns shows 1 if power supply is in CC mode

property cv_mode: hvl_ccb.dev.fug.FuGDigitalVal
Returns shows 1 if power supply is in CV mode

property digital_control: hvl_ccb.dev.fug.FuGDigitalVal
Returns shows 1 if power supply is digitally controlled

property on: hvl_ccb.dev.fug.FuGDigitalVal
Returns shows 1 if power supply ON

property reg_3: hvl_ccb.dev.fug.FuGDigitalVal
For special applications.

Returns input from bit 3-REG

property x_stat: hvl_ccb.dev.fug.FuGPolarities
Returns polarity of HVPS with polarity reversal

class FuGProbusVDORegisters(fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups)
Bases: object

Digital outputs acc. 4.2.2

property out: Union[int, hvl_ccb.dev.fug.FuGDigitalVal]
Status of the output according to the last setting. This can differ from the actual state if output should only pulse.

Returns FuGDigitalVal

property status: `hvl_ccb.dev.fug.FuGDigitalVal`

Returns the actual value of output. This can differ from the set value if pulse function is used.

Returns FuGDigitalVal

class FuGProbusVMonitorRegisters(*fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Analog monitors acc. 4.2.3

property adc_mode: `hvl_ccb.dev.fug.FuGMonitorModes`

The programmed resolution and integration time of the AD converter

Returns FuGMonitorModes

property value: `float`

Value from the monitor.

Returns a float value in V or A

property value_raw: `float`

uncalibrated raw value from AD converter

Returns float value from ADC

class FuGProbusVRegisterGroups(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)

Bases: `hvl_ccb.utils.enum.NameEnum`

An enumeration.

CONFIG = 'K'

INPUT = 'D'

MONITOR_I = 'M1'

MONITOR_V = 'M0'

OUTPUTONCMD = 'BON'

OUTPUTX0 = 'B0'

OUTPUTX1 = 'B1'

OUTPUTX2 = 'B2'

OUTPUTXCMD = 'BX'

SETCURRENT = 'S1'

SETVOLTAGE = 'S0'

class FuGProbusVSetRegisters(*fug, super_register: hvl_ccb.dev.fug.FuGProbusVRegisterGroups*)

Bases: object

Setvalue control acc. 4.2.1 for the voltage and the current output

property actualsetvalue: `float`

The actual valid set value, which depends on the ramp function.

Returns actual valid set value

property high_resolution: `hvl_ccb.dev.fug.FuGDigitalVal`

Status of the high resolution mode of the output.

Return 0 normal operation

Return 1 High Res. Mode

property rampmode: *hvl_ccb.dev.fug.FuGRampModes*

The set ramp mode to control the setvalue.

Returns the mode of the ramp as instance of FuGRampModes

property ramprate: *float*

The set ramp rate in V/s.

Returns ramp rate in V/s

property rampstate: *hvl_ccb.dev.fug.FuGDigitalVal*

Status of ramp function.

Return 0 if final setvalue is reached

Return 1 if still ramping up

property setvalue: *float*

For the voltage or current output this setvalue was programmed.

Returns the programmed setvalue

class FuGRampModes(*value*)

Bases: enum.IntEnum

An enumeration.

FOLLOWRAMP = 1

Follow the ramp up- and downwards

IMMEDIATELY = 0

Standard mode: no ramp

ONLYUPWARDSOFFTOZERO = 4

Follow the ramp up- and downwards, if output is OFF set value is zero

RAMPUPWARDS = 2

Follow the ramp only upwards, downwards immediately

SPECIALRAMPUPWARDS = 3

Follow a special ramp function only upwards

class FuGReadbackChannels(*value*)

Bases: enum.IntEnum

An enumeration.

CURRENT = 1

FIRMWARE = 5

RATEDCURRENT = 4

RATEDVOLTAGE = 3

SN = 6

STATUSBYTE = 2

VOLTAGE = 0

class FuGSerialCommunication(*configuration*)

Bases: *hvl_ccb.comm.serial.SerialCommunication*

Specific communication protocol implementation for FuG power supplies. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(command: str) → str

Send a command to the interface and handle the status message. Eventually raises an exception.

Parameters **command** – Command to send

Raises **FuGError** – if the connection is broken or the error from the power source itself

Returns Answer from the interface or empty string

```
class FuGSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
                                    encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty:
                                    Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int =
                                    10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
                                    <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,
                                    hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                    <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,
                                    hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int,
                                    float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for FuG power supplies is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

default_n_attempts_read_text_nonempty: int = 10

default number of attempts to read a non-empty text

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

FuG does not use parity

```
classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
    FuG uses one stop bit

terminator: bytes = b'\n'
    The terminator is LF

timeout: Union[int, float] = 3
    use 3 seconds timeout as default

wait_sec_read_text_nonempty: Union[int, float] = 0.5
    default time to wait between attempts of reading a non-empty text

class FuGTerminators(value)
Bases: enum.IntEnum

An enumeration.


CR = 3

CRLF = 0

LF = 2

LFCR = 1
```

hvl_ccb.dev.heinzinger

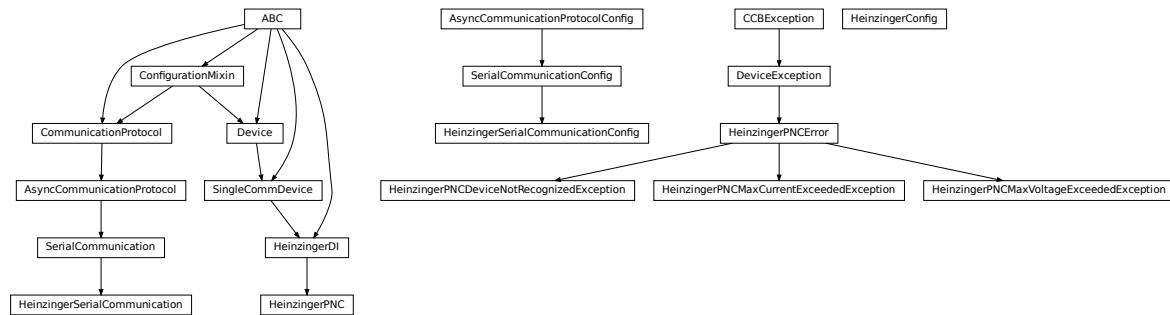

Device classes for Heinzinger Digital Interface I/II and Heinzinger PNC power supply.

The Heinzinger Digital Interface I/II is used for many Heinzinger power units. Manufacturer homepage: <https://www.heinzinger.com/products/accessories-and-more/digital-interfaces/>

The Heinzinger PNC series is a series of high voltage direct current power supplies. The class HeinzingerPNC is tested with two PNChp 60000-1neg and a PNChp 1500-1neg. Check the code carefully before using it with other PNC devices, especially PNC3p or PNCCap. Manufacturer homepage: <https://www.heinzinger.com/products/high-voltage/universal-high-voltage-power-supplies/>

```
class HeinzingerConfig(default_number_of_recordings: Union[int,
    hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum] = 1,
    number_of_decimals: int = 6, wait_sec_stop_commands: Union[int, float] = 0.5)
Bases: object
```

Device configuration dataclass for Heinzinger power supplies.

```
class RecordingsEnum(value)
    Bases: enum.IntEnum

    An enumeration.

    EIGHT = 8
    FOUR = 4
    ONE = 1
    SIXTEEN = 16
    TWO = 2

clean_values()

default_number_of_recordings: Union[int,
    hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum] = 1

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.
```

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

number_of_decimals: int = 6

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

wait_sec_stop_commands: Union[int, float] = 0.5

Time to wait after subsequent commands during stop (in seconds)

class HeinzingerDI(*com*, *dev_config=None*)

Bases: *hvl_ccb.dev.base.SingleCommDevice*, abc.ABC

Heinzinger Digital Interface I/II device class

Sends basic SCPI commands and reads the answer. Only the standard instruction set from the manual is implemented.

```
class OutputStatus(value)
    Bases: enum.IntEnum

    Status of the voltage output

    OFF = 0
    ON = 1
    UNKNOWN = -1

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

get_current() → float
    Queries the set current of the Heinzinger PNC (not the measured current!).

    Raises SerialCommunicationIOError – when communication port is not opened

get_interface_version() → str
    Queries the version number of the digital interface.

    Raises SerialCommunicationIOError – when communication port is not opened

get_number_of_recordings() → int
    Queries the number of recordings the device is using for average value calculation.

    Returns int number of recordings

    Raises SerialCommunicationIOError – when communication port is not opened

get_serial_number() → str
    Ask the device for its serial number and returns the answer as a string.

    Returns string containing the device serial number

    Raises SerialCommunicationIOError – when communication port is not opened

get_voltage() → float
    Queries the set voltage of the Heinzinger PNC (not the measured voltage!).

    Raises SerialCommunicationIOError – when communication port is not opened

measure_current() → float
    Ask the Device to measure its output current and return the measurement result.

    Returns measured current as float

    Raises SerialCommunicationIOError – when communication port is not opened

measure_voltage() → float
    Ask the Device to measure its output voltage and return the measurement result.

    Returns measured voltage as float

    Raises SerialCommunicationIOError – when communication port is not opened

output_off() → None
    Switch DC voltage output off and updates the output status.

    Raises SerialCommunicationIOError – when communication port is not opened
```

output_on() → None
 Switch DC voltage output on and updates the output status.

Raises `SerialCommunicationIOError` – when communication port is not opened

property output_status: `hvl_ccb.dev.heinzinger.HeinzingerDI.OutputStatus`

reset_interface() → None
 Reset of the digital interface; only Digital Interface I: Power supply is switched to the Local-Mode (Manual operation)

Raises `SerialCommunicationIOError` – when communication port is not opened

set_current(value: Union[int, float]) → None
 Sets the output current of the Heinzinger PNC to the given value.

Parameters `value` – current expressed in `self.unit_current`

Raises `SerialCommunicationIOError` – when communication port is not opened

set_number_of_recordings(value: Union[int, hvl_ccb.dev.heinzinger.HeinzingerConfig.RecordingsEnum]) → None
 Sets the number of recordings the device is using for average value calculation. The possible values are 1, 2, 4, 8 and 16.

Raises `SerialCommunicationIOError` – when communication port is not opened

set_voltage(value: Union[int, float]) → None
 Sets the output voltage of the Heinzinger PNC to the given value.

Parameters `value` – voltage expressed in `self.unit_voltage`

Raises `SerialCommunicationIOError` – when communication port is not opened

abstract start()
 Opens the communication protocol.

Raises `SerialCommunicationIOError` – when communication port cannot be opened.

stop() → None
 Stop the device. Closes also the communication protocol.

class HeinzingerPNC(com, dev_config=None)
 Bases: `hvl_ccb.dev.heinzinger.HeinzingerDI`

Heinzinger PNC power supply device class.

The power supply is controlled over a Heinzinger Digital Interface I/II

class UnitCurrent(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
 Bases: `hvl_ccb.utils.enum.AutoNumberNameEnum`

An enumeration.

A = 3

UNKNOWN = 1

mA = 2

class UnitVoltage(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
 Bases: `hvl_ccb.utils.enum.AutoNumberNameEnum`

An enumeration.

UNKNOWN = 1

```
V = 2
kV = 3

identify_device() → None
    Identify the device nominal voltage and current based on its serial number.

    Raises SerialCommunicationIOError – when communication port is not opened

property max_current: Union[int, float]
property max_current_hardware: Union[int, float]
property max_voltage: Union[int, float]
property max_voltage_hardware: Union[int, float]

set_current(value: Union[int, float]) → None
    Sets the output current of the Heinzinger PNC to the given value.

    Parameters value – current expressed in self.unit_current

    Raises SerialCommunicationIOError – when communication port is not opened

set_voltage(value: Union[int, float]) → None
    Sets the output voltage of the Heinzinger PNC to the given value.

    Parameters value – voltage expressed in self.unit_voltage

    Raises SerialCommunicationIOError – when communication port is not opened

start() → None
    Opens the communication protocol and configures the device.

property unit_current: hvl_ccb.dev.heinzinger.HeinzingerPNC.UnitCurrent
property unit_voltage: hvl_ccb.dev.heinzinger.HeinzingerPNC.UnitVoltage

exception HeinzingerPNCDriverNotRecognizedException
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

    Error indicating that the serial number of the device is not recognized.

exception HeinzingerPNCError
    Bases: hvl_ccb.dev.base.DeviceException

    General error with the Heinzinger PNC voltage source.

exception HeinzingerPNCMaxCurrentExceededError
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

    Error indicating that program attempted to set the current to a value exceeding ‘max_current’.

exception HeinzingerPNCMaxVoltageExceededError
    Bases: hvl_ccb.dev.heinzinger.HeinzingerPNCError

    Error indicating that program attempted to set the voltage to a value exceeding ‘max_voltage’.

class HeinzingerSerialCommunication(configuration)
    Bases: hvl_ccb.comm.serial.SerialCommunication

    Specific communication protocol implementation for Heinzinger power supplies. Already predefines device-specific protocol parameters in config.

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class
```

```
class HeinzingerSerialCommunicationConfig(terminator: bytes = b'\n', encoding: str = 'utf-8',
    encoding_error_handling: str = 'strict',
    wait_sec_read_text_nonempty: Union[int, float] = 0.5,
    default_n_attempts_read_text_nonempty: int = 40, port:
    Union[str, NoneType] = None, baudrate: int = 9600, parity:
    Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] =
    <SerialCommunicationParity.NONE: 'N'>, stopbits:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationStopbits] =
    <SerialCommunicationStopbits.ONE: 1>, bytesize:
    Union[int,
    hvl_ccb.comm.serial.SerialCommunicationBytesize] =
    <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
    Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for Heinzinger power supplies is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

default_n_attempts_read_text_nonempty: int = 40

increased to 40 default number of attempts to read a non-empty text

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

Heinzinger does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1

Heinzinger uses one stop bit


terminator: bytes = b'\n'

The terminator is LF

```
timeout: Union[int, float] = 3
use 3 seconds timeout as default

wait_sec_read_text_nonempty: Union[int, float] = 0.5
default time to wait between attempts of reading a non-empty text
```

hvl_ccb.dev.labjack

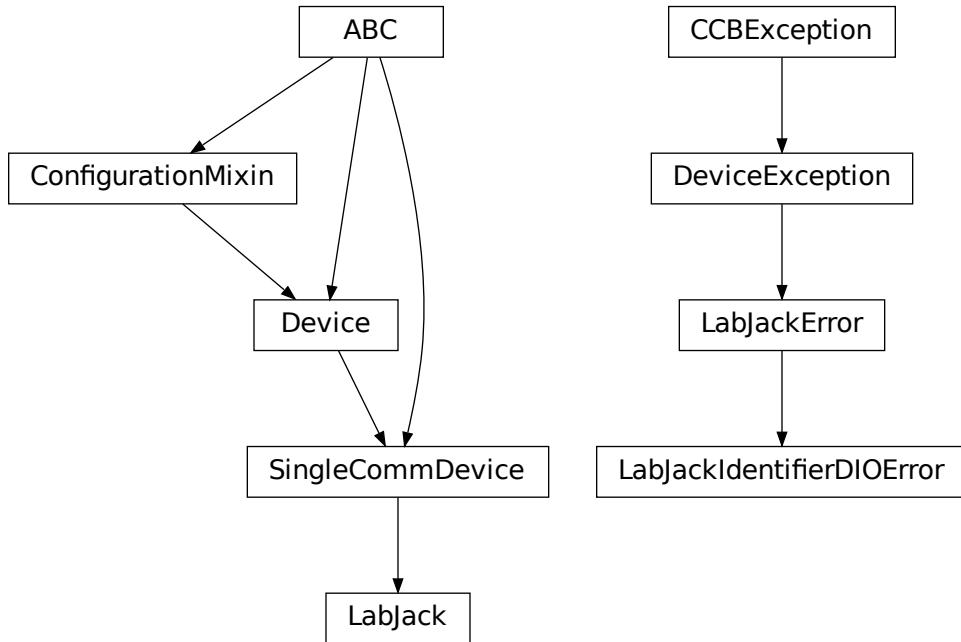

A
LabJack T-series devices wrapper around the LabJack's LJM Library; see <https://labjack.com/ljm> . The wrapper was originally developed and tested for a LabJack T7-PRO device.

Extra installation

To use this LabJack T-series devices wrapper:

1. install the hvl_ccb package with a labjack extra feature:

```
$ pip install "hvl_ccb[labjack]"
```

this will install the Python bindings for the library.

2. install the library - follow instruction in <https://labjack.com/support/software/installers/ljm> .

```
class LabJack(com, dev_config=None)
    Bases: hvl_ccb.dev.base.SingleCommDevice

    LabJack Device.
```

This class is tested with a LabJack T7-Pro and should also work with T4 and T7 devices communicating through the LJM Library. Other or older hardware versions and variants of LabJack devices are not supported.

```
class AInRange(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl\_ccb.utils.enum.StrEnumBase
```

An enumeration.

ONE = 1.0

ONE_HUNDREDTH = 0.01

ONE_TENTH = 0.1

TEN = 10.0

property value: float

```
class BitLimit(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: aenum.IntEnum
```

Maximum integer values for clock settings

THIRTY_TWO_BIT = 4294967295

```
class CalMicroAmpere(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
    Bases: aenum.Enum
```

Pre-defined microampere (uA) values for calibration current source query.

TEN = '10uA'

TWO_HUNDRED = '200uA'

```
class CjcType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl\_ccb.utils.enum.NameEnum
```

CJC slope and offset

internal = (1, 0)

lm34 = (55.56, 255.37)

```
class ClockFrequency(value=<no_arg>, names=None, module=None, type=None, start=1,
                           boundary=None)
    Bases: aenum.IntEnum
```

Available clock frequencies, in Hz

FIVE_MHZ = 5000000

FORTY_MHZ = 40000000

MAXIMUM = 80000000

MINIMUM = 312500

TEN_MHZ = 10000000

TWELVE_HUNDRED_FIFTY_KHZ = 1250000

TWENTY_FIVE_HUNDRED_KHZ = 2500000

TWENTY_MHZ = 20000000

DIOChannel

alias of `hvl_ccb._dev.labjack.TSeriesDIOChannel`

```
class DIOStatus(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: aenum.IntEnum
State of a digital I/O channel.

HIGH = 1
LOW = 0

class DeviceType(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
Bases: hvl_ccb.utils.enum.AutoNumberNameEnum
LabJack device types.

Can be also looked up by ambiguous Product ID (p_id) or by instance name: `python
LabJackDeviceType(4) is LabJackDeviceType('T4')` 

ANY = 1
T4 = 2
T7 = 3
T7_PRO = 4

@classmethod get_by_p_id(p_id: int) → Union[hvl_ccb._dev.labjack.DeviceType,
List[hvl_ccb._dev.labjack.DeviceType]]
Get LabJack device type instance via LabJack product ID.

Note: Product ID is not unambiguous for LabJack devices.
Parameters p_id – Product ID of a LabJack device
Returns Instance or list of instances of LabJackDeviceType
Raises ValueError – when Product ID is unknown

class TemperatureUnit(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
Temperature unit (to be returned)

C = 1
F = 2
K = 0

class ThermocoupleType(value=<no_arg>, names=None, module=None, type=None, start=1,
boundary=None)
Bases: hvl_ccb.utils.enum.NameEnum
Thermocouple type; NONE means disable thermocouple mode.

C = 30
E = 20
J = 21
K = 22
NONE = 0
PT100 = 40
PT1000 = 42
PT500 = 41
```

R = 23**S = 25****T = 24**

config_high_pulse(*address*: Union[str, *hvl_ccb._dev.labjack.TSeriesDIOChannel*], *t_start*: Union[int, float], *t_width*: Union[int, float], *n_pulses*: int = 1) → None

Configures one FIO channel to send a timed HIGH pulse. Configure multiple channels to send pulses with relative timing accuracy. Times have a maximum resolution of 1e-7 seconds @ 10 MHz. :param address: FIO channel: [T7] FIO0;2;3;4;5. [T4] FIO6;7. :raises LabJackError if address is not supported. :param t_start: pulse start time in seconds. :raises ValueError: if t_start is negative or would exceed the clock period. :param t_width: duration of high pulse, in seconds. :raises ValueError: if t_width is negative or would exceed the clock period. :param n_pulses: number of pulses to be sent; single pulse default. :raises TypeError if n_pulses is not of type int. :raises ValueError if n_pulses is negative or exceeds the 32bit limit.

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

disable_pulses(**addresses*: Optional[Union[str, *hvl_ccb._dev.labjack.TSeriesDIOChannel*]]) → None

Disable previously configured pulse channels. :param addresses: tuple of FIO addresses. All channels disabled if no argument is passed.

enable_clock(*clock_enabled*: bool) → None

Enable/disable LabJack clock to configure or send pulses. :param clock_enabled: True -> enable, False -> disable. :raises TypeError: if clock_enabled is not of type bool

get_ain(**channels*: int) → Union[float, Sequence[float]]

Read currently measured value (voltage, resistance, ...) from one or more of analog inputs.

Parameters **channels** – AIN number or numbers (0..254)

Returns the read value (voltage, resistance, ...) as float or tuple of them in case multiple channels given

get_cal_current_source(*name*: Union[str, *CalMicroAmpere*]) → float

This function will return the calibration of the chosen current source, this ist not a measurement!

The value was stored during fabrication.

Parameters **name** – ‘200uA’ or ‘10uA’ current source

Returns calibration of the chosen current source in ampere

get_clock() → Dict[str, Union[int, float]]

Return clock settings read from LabJack.

get_digital_input(*address*: Union[str, *hvl_ccb._dev.labjack.TSeriesDIOChannel*]) → *hvl_ccb.dev.labjack.LabJack.DIOSStatus*

Get the value of a digital input.

allowed names for T7 (Pro): FIO0 - FIO7, EIO0 - EIO 7, CIO0- CIO3, MIO0 - MIO2 :param address: name of the output -> ‘FIO0’ :return: HIGH when *address* DIO is high, and LOW when *address* DIO is low

get_product_id() → int

This function returns the product ID reported by the connected device.

Attention: returns 7 for both T7 and T7-Pro devices!

Returns integer product ID of the device

get_product_name(*force_query_id=False*) → str
This function will return the product name based on product ID reported by the device.
Attention: returns “T7” for both T7 and T7-Pro devices!

Parameters **force_query_id** – boolean flag to force *get_product_id* query to device instead of using cached device type from previous queries.

Returns device name string, compatible with *LabJack.DeviceType*

get_product_type(*force_query_id: bool = False*) → *hvl_ccb._dev.labjack.DeviceType*
This function will return the device type based on reported device type and in case of unambiguity based on configuration of device’s communication protocol (e.g. for “T7” and “T7_PRO” devices), or, if not available first matching.

Parameters **force_query_id** – boolean flag to force *get_product_id* query to device instead of using cached device type from previous queries.

Returns *DeviceType* instance

Raises *LabJackIdentifierDIOError* – when read Product ID is unknown

get_sbus_rh(*number: int*) → float
Read the relative humidity value from a serial SBUS sensor.

Parameters **number** – port number (0..22)

Returns relative humidity in %RH

get_sbus_temp(*number: int*) → float
Read the temperature value from a serial SBUS sensor.

Parameters **number** – port number (0..22)

Returns temperature in Kelvin

get_serial_number() → int
Returns the serial number of the connected LabJack.

Returns Serial number.

read_resistance(*channel: int*) → float
Read resistance from specified channel.

Parameters **channel** – channel with resistor

Returns resistance value with 2 decimal places

read_thermocouple(*pos_channel: int*) → float
Read the temperature of a connected thermocouple.

Parameters **pos_channel** – is the AIN number of the positive pin

Returns temperature in specified unit

send_pulses(**addresses: Union[str, hvl_ccb._dev.labjack.TSeriesDIOChannel]*) → None
Sends pre-configured pulses for specified addresses. :param *addresses*: tuple of FIO addresses :raises LabJackError if an address has not been configured.

set_ain_differential(*pos_channel: int, differential: bool*) → None
Sets an analog input to differential mode or not. T7-specific: For base differential channels, positive must be even channel from 0-12 and negative must be positive+1. For extended channels 16-127, see Mux80 datasheet.

Parameters

- **pos_channel** – is the AIN number (0..12)
- **differential** – True or False

Raises `LabJackError` – if parameters are unsupported

set_ain_range(*channel*: int, *vrange*: Union[Real, AInRange]) → None

Set the range of an analog input port.

Parameters

- **channel** – is the AIN number (0..254)
- **vrange** – is the voltage range to be set

set_ain_resistance(*channel*: int, *vrange*: Union[Real, AInRange], *resolution*: int) → None

Set the specified channel to resistance mode. It utilized the 200uA current source of the LabJack.

Parameters

- **channel** – channel that should measure the resistance
- **vrange** – voltage range of the channel
- **resolution** – resolution index of the channel T4: 0-5, T7: 0-8, T7-Pro 0-12

set_ain_resolution(*channel*: int, *resolution*: int) → None

Set the resolution index of an analog input port.

Parameters

- **channel** – is the AIN number (0..254)
- **resolution** – is the resolution index within 0...`get_product_type().ain_max_resolution` range; 0 will set the resolution index to default value.

set_ain_thermocouple(*pos_channel*: int, *thermocouple*: Union[None, str, ThermocoupleType],
cjc_address: int = 60050, *cjc_type*: Union[str, CjcType] = CjcType.internal,
vrange: Union[Real, AInRange] = AInRange.ONE_HUNDREDTH, *resolution*: int
= 10, *unit*: Union[str, TemperatureUnit] = TemperatureUnit.K) → None

Set the analog input channel to thermocouple mode.

Parameters

- **pos_channel** – is the analog input channel of the positive part of the differential pair
- **thermocouple** – None to disable thermocouple mode, or string specifying the thermocouple type
- **cjc_address** – modbus register address to read the CJC temperature
- **cjc_type** – determines cjc slope and offset, ‘internal’ or ‘lm34’
- **vrange** – measurement voltage range
- **resolution** – resolution index (T7-Pro: 0-12)
- **unit** – is the temperature unit to be returned (‘K’, ‘C’ or ‘F’)

Raises `LabJackError` – if parameters are unsupported

set_analog_output(*channel*: int, *value*: Union[int, float]) → None

Set the voltage of a analog output port

Parameters

- **channel** – DAC channel number 1/0

- **value** – The output voltage value 0-5 Volts int/float

set_clock(*clock_frequency*: Union[Number, ClockFrequency] = 10000000, *clock_period*: Number = 1) → None

Configure LabJack clock for pulse out feature. :param *clock_frequency*: clock frequency in Hz; default 10 MHz for base 10. :raises ValueError: if *clock_frequency* is not allowed (see ClockFrequency). :param *clock_period*: clock roll time in seconds; default 1s, 0 for max. :raises ValueError: if *clock_period* exceeds the 32bit tick limit. Clock period determines pulse spacing when using multi-pulse settings. Ensure period exceeds maximum intended pulse end time.

set_digital_output(*address*: str, *state*: Union[int, DIOStatus]) → None

Set the value of a digital output.

Parameters

- **address** – name of the output -> ‘FIO0’
- **state** – state of the output -> *DIOStatus* instance or corresponding *int* value

start() → None

Start the Device.

stop() → None

Stop the Device.

exception LabJackError

Bases: *hvl_ccb.dev.base.DeviceException*


General Exception for the LabJack device.

exception LabJackIdentifierDIOError

Bases: *hvl_ccb.dev.labjack.LabJackError*

Error indicating a wrong DIO identifier

hvl_ccb.dev.lauda

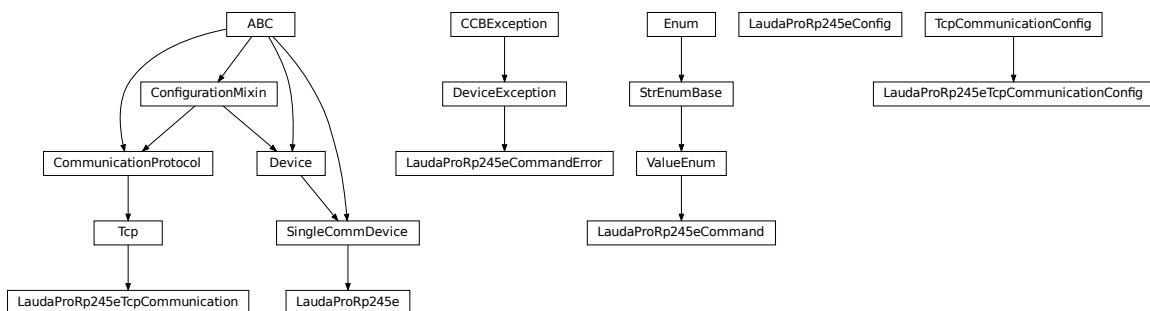

Device class for controlling a Lauda PRO RP245E, circulation chiller over TCP.

class LaudaProRp245e(*com*, *dev_config*=None)

Bases: *hvl_ccb.dev.base.SingleCommDevice*

Lauda RP245E circulation chiller class.

static config_cls() → Type[*hvl_ccb.dev.lauda.LaudaProRp245eConfig*]

Return the default configdataclass class.

Returns a reference to the default configdataclass class

continue_ramp() → str
Continue current ramp program.

Returns reply of the device to the last call of “query”

static default_com_cls() → Type[hvl_ccb.dev.lauda.LaudaProRp245eTcpCommunication]
Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

get_bath_temp() → float
:return : float value of measured lauda bath temp in °C

get_device_type() → str
:return : Connected Lauda device type (for connection/com test)

pause() → str
Stop temperature control and pump.

Returns reply of the device to the last call of “query”

pause_ramp() → str
Pause current ramp program.

Returns reply of the device to the last call of “query”

reset_ramp() → str
Delete all segments from current ramp program.

Returns reply of the device to the last call of “query”

run() → str
Start temperature control & pump.

Returns reply of the device to the last call of “query”

set_control_mode(mod: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = ExtControlModeEnum.INTERNAL) → str
Define control mode. 0 = INTERNAL (control bath temp), 1 = EXPT100 (pt100 attached to chiller), 2 = ANALOG, 3 = SERIAL, 4 = USB, 5 = ETH (to be used when passing the ext. temp. via ethernet) (temperature then needs to be passed every second, when not using options 3, 4, or 5)

Parameters **mod** – temp control mode (control internal temp or external temp).

Returns reply of the device to the last call of “query” (“OK”, if command was recognized”)

set_external_temp(external_temp: float = 20.0) → str
Pass value of external controlled temperature. Should be done every second, when control of external temperature is active. Has to be done right before control of external temperature is activated.

Parameters **external_temp** – current value of external temperature to be controlled.

Returns reply of the device to the last call of “query”

set_pump_level(pump_level: int = 6) → str
Set pump level Raises ValueError, if pump level is invalid.

Parameters **pump_level** – pump level.

Returns reply of the device to the last call of “query”

set_ramp_iterations(num: int = 1) → str
Define number of ramp program cycles.

Parameters `num` – number of program cycles to be performed.

Returns reply of the device to the last call of “query”

set_ramp_program(`program: int = 1`) → str

Define ramp program for following ramp commands. Raises ValueError if maximum number of ramp programs (5) is exceeded.

Parameters `program` – Number of ramp program to be activated for following commands.

Returns reply of the device to the last call of “query”

set_ramp_segment(`temp: float = 20.0, dur: int = 0, tol: float = 0.0, pump: int = 6`) → str

Define segment of current ramp program - will be attached to current program. Raises ValueError, if pump level is invalid.

Parameters

- `temp` – target temperature of current ramp segment
- `dur` – duration in minutes, in which target temperature should be reached
- `tol` – tolerance at which target temperature should be reached (for 0.00, next segment is started after dur has passed).
- `pump` – pump level to be used for this program segment.

Returns reply of the device to the last call of “query”

set_temp_set_point(`temp_set_point: float = 20.0`) → str

Define temperature set point

Parameters `temp_set_point` – temperature set point.

Returns reply of the device to the last call of “query”

start() → None

Start this device.

start_ramp() → str

Start current ramp program.

Returns reply of the device to the last call of “query”

stop() → None

Stop this device. Disables access and closes the communication protocol.

stop_ramp() → str

Stop current ramp program.

Returns reply of the device to the last call of “query”

validate_pump_level(`level: int`)

Validates pump level. Raises ValueError, if pump level is incorrect. :param level: pump level, integer

class LaudaProRp245eCommand(`value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None`)

Bases: `hvl_ccb.utils.enum.ValueEnum`

Commands for Lauda PRO RP245E Chiller Command strings most often need to be complimented with a parameter (attached as a string) before being sent to the device. Commands implemented as defined in “Lauda Betriebsanleitung fuer PRO Badthermostate und Umwaelzthermostate” pages 42 - 49

BATH_TEMP = 'IN_PV_00'

Request internal bath temperature

```

COM_TIME_OUT = 'OUT_SP_08_'
    Define communication time out

CONT_MODE = 'OUT_MODE_01_'
    Set control mode 1=internal, 2=ext. analog, 3=ext. serial, 4=USB, 5=ethernet

DEVICE_TYPE = 'TYPE'
    Request device type

EXTERNAL_TEMP = 'OUT_PV_05_'
    Pass on external controlled temperature

LOWER_TEMP = 'OUT_SP_05_'
    Define lower temp limit

OPERATION_MODE = 'OUT_SP_02_'
    Define operation mode

PUMP_LEVEL = 'OUT_SP_01_'
    Define pump level 1-8

RAMP_CONTINUE = 'RMP_CONT'
    Continue a paused ramp program

RAMP_DELETE = 'RMP_RESET'
    Reset a selected ramp program

RAMP_ITERATIONS = 'RMP_OUT_02_'
    Define how often a ramp program should be iterated

RAMP_PAUSE = 'RMP_PAUSE'
    Pause a selected ramp program

RAMP_SELECT = 'RMP_SELECT_'
    Select a ramp program (target for all further ramp commands)

RAMP_SET = 'RMP_OUT_00_'
    Define parameters of a selected ramp program

RAMP_START = 'RMP_START'
    Start a selected ramp program

RAMP_STOP = 'RMP_STOP'
    Stop a running ramp program

START = 'START'
    Start temp control (pump and heating/cooling)

STOP = 'STOP'
    Stop temp control (pump and heating/cooling)

TEMP_SET_POINT = 'OUT_SP_00_'
    Define temperature set point

UPPER_TEMP = 'OUT_SP_04_'
    Define upper temp limit

build_str(param: str = "", terminator: str = '\r\n')
    Build a command string for sending to the device

```

Parameters

- **param** – Command's parameter given as string
- **terminator** – Command's terminator

Returns Command's string with a parameter and terminator

exception LaudaProRp245eCommandError

Bases: `hvl_ccb.dev.base.DeviceException`

Exception raised when an error is returned upon a command.

class LaudaProRp245eConfig(`temp_set_point_init: Union[int, float] = 20.0, pump_init: int = 6, upper_temp: Union[int, float] = 80.0, lower_temp: Union[int, float] = -55.0, com_time_out: Union[int, float] = 0, max_pump_level: int = 8, max_pr_number: int = 5, operation_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.OperationModeEnum] = OperationModeEnum.AUTO, control_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = ExtControlModeEnum.INTERNAL)`)

Bases: `object`

Configuration for the Lauda RP245E circulation chiller.

class ExtControlModeEnum(`value`)

Bases: `enum.IntEnum`

Source for definition of external, controlled temperature (option 2, 3 and 4 are not available with current configuration of the Lauda RP245E, add-on hardware would required)

`ANALOG = 2`

`ETH = 5`

`EXPT100 = 1`

`INTERNAL = 0`

`SERIAL = 3`

`USB = 4`

class OperationModeEnum(`value`)

Bases: `enum.IntEnum`

Operation Mode (Cooling OFF/Cooling On/AUTO - set to AUTO)

`AUTO = 2`

Automatically select heating/cooling

`COOLOFF = 0`

`COOLON = 1`

`clean_values() → None`

`com_time_out: Union[int, float] = 0`

Communication time out (0 = OFF)

`control_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.ExtControlModeEnum] = 0`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- `fieldname` – name of the field

- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
 Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

lower_temp: Union[int, float] = -55.0
 Lower temperature limit (safe for Galden HT135 cooling liquid)

max_pr_number: int = 5
 Maximum number of ramp programs that can be stored in the memory of the chiller

max_pump_level: int = 8
 Highest pump level of the chiller

operation_mode: Union[int, hvl_ccb.dev.lauda.LaudaProRp245eConfig.OperationModeEnum] = 2

classmethod optional_defaults() → Dict[str, object]
 Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

pump_init: int = 6
 Default pump Level

classmethod required_keys() → Sequence[str]
 Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

temp_set_point_init: Union[int, float] = 20.0
 Default temperature set point

upper_temp: Union[int, float] = 80.0
 Upper temperature limit (safe for Galden HT135 cooling liquid)

class LaudaProRp245eTcpCommunication(configuration)
 Bases: *hvl_ccb.comm.tcp.Tcp*

Implements the Communication Protocol for Lauda PRO RP245E TCP connection.

close() → None
 Close the Lauda PRO RP245E TCP connection.

static config_cls() → Type[hvl_ccb.dev.lauda.LaudaProRp245eTcpCommunicationConfig]
 Return the default configdataclass class.

Returns a reference to the default configdataclass class

open() → None
 Open the Lauda PRO RP245E TCP connection.

Raises *LaudaProRp245eCommandError* – if the connection fails.

query_command(command: hvl_ccb.dev.lauda.LaudaProRp245eCommand, param: str = "") → str
 Send and receive function. E.g. to be used when setting/changing device setting. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None

read() → str

Receive value function. :return: reply from device as a string, the terminator, as well as the ‘OK’ stripped from the reply to make it directly useful as a value (e.g. in case the internal bath temperature is requested)

write_command(command: hvl_ccb.dev.lauda.LaudaProRp245eCommand, param: str = "") → None

Send command function. :param command: first part of command string, defined in *LaudaProRp245eCommand* :param param: second part of command string, parameter (by default '') :return: None

class LaudaProRp245eTcpCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], port: int = 54321, bufsize: int = 1024, wait_sec_pre_read_or_write: Union[int, float] = 0.005, terminator: str = '\r\n')

Bases: *hvl_ccb.comm.tcp.TcpCommunicationConfig*

Configuration dataclass for *LaudaProRp245eTcpCommunication*.

clean_values() → None

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.


Returns a list of strings containing all required keys.

terminator: str = '\r\n'

The terminator character

wait_sec_pre_read_or_write: Union[int, float] = 0.005

Delay time between commands in seconds

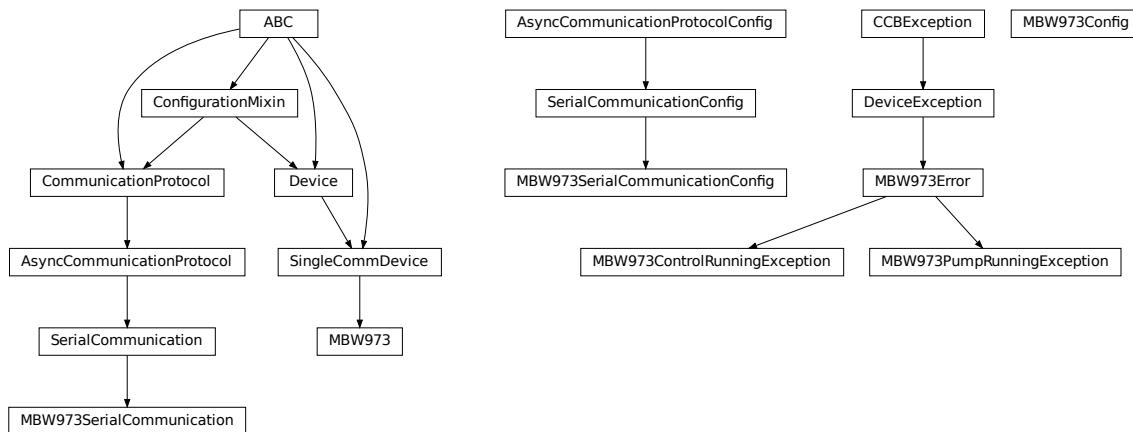hvl_ccb.dev.mbw973

Device class for controlling a MBW 973 SF6 Analyzer over a serial connection.

The MBW 973 is a gas analyzer designed for gas insulated switchgear and measures humidity, SF6 purity and SO2 contamination in one go. Manufacturer homepage: <https://www.mbw.ch/products/sf6-gas-analysis/973-sf6-analyzer/>

```

class MBW973(com, dev_config=None)
  Bases: hvl_ccb.dev.base.SingleCommDevice

  MBW 973 dew point mirror device class.

  static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

  static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

  is_done() → bool
    Poll status of the dew point mirror and return True, if all measurements are done.

    Returns True, if all measurements are done; False otherwise.

  Raises SerialCommunicationIOError – when communication port is not opened

  read(cast_type: Type = <class 'str'>)
    Read value from self.com and cast to cast_type. Raises ValueError if read text (str) is not convertible to cast_type, e.g. to float or to int.

    Returns Read value of cast_type type.

  read_float() → float
    Convenience wrapper for self.read(), with typing hint for return value.

    Returns Read float value.

  read_int() → int
    Convenience wrapper for self.read(), with typing hint for return value.
  
```

Returns Read *int* value.

read_measurements() → Dict[str, float]

Read out measurement values and return them as a dictionary.

Returns Dictionary with values.

Raises *SerialCommunicationIOError* – when communication port is not opened

set_measuring_options(humidity: bool = True, sf6_purity: bool = False) → None

Send measuring options to the dew point mirror.

Parameters

- **humidity** – Perform humidity test or not?

- **sf6_purity** – Perform SF6 purity test or not?

Raises *SerialCommunicationIOError* – when communication port is not opened

start() → None

Start this device. Opens the communication protocol and retrieves the set measurement options from the device.

Raises *SerialCommunicationIOError* – when communication port cannot be opened.

start_control() → None

Start dew point control to acquire a new value set.

Raises *SerialCommunicationIOError* – when communication port is not opened

stop() → None

Stop the device. Closes also the communication protocol.

write(value) → None

Send *value* to *self.com*.

Parameters **value** – Value to send, converted to *str*.

Raises *SerialCommunicationIOError* – when communication port is not opened

class MBW973Config(polling_interval: Union[int, float] = 2)

Bases: *object*

Device configuration dataclass for MBW973.

clean_values()

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field

- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

polling_interval: Union[int, float] = 2

Polling period for *is_done* status queries [in seconds].

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception MBW973ControlRunningException

Bases: *hvl_ccb.dev.mbw973.MBW973Error*

Error indicating there is still a measurement running, and a new one cannot be started.

exception MBW973Error

Bases: *hvl_ccb.dev.base.DeviceException*

General error with the MBW973 dew point mirror device.

exception MBW973PumpRunningException

Bases: *hvl_ccb.dev.mbw973.MBW973Error*

Error indicating the pump of the dew point mirror is still recovering gas, unable to start a new measurement.

class MBW973SerialCommunication(configuration)

Bases: *hvl_ccb.comm.serial.SerialCommunication*

Specific communication protocol implementation for the MBW973 dew point mirror. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class MBW973SerialCommunicationConfig(terminator: bytes = b'\r', encoding: str = 'utf-8',

encoding_error_handling: str = 'strict',

wait_sec_read_text_nonempty: Union[int, float] = 0.5,

default_n_attempts_read_text_nonempty: int = 10, port: Union[str,

NoneType] = None, baudrate: int = 9600, parity: Union[str,

hvl_ccb.comm.serial.SerialCommunicationParity] =

<SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int,

hvl_ccb.comm.serial.SerialCommunicationStopbits] =

<SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int,

hvl_ccb.comm.serial.SerialCommunicationBytesize] =

<SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:

Union[int, float] = 3)

Bases: *hvl_ccb.comm.serial.SerialCommunicationConfig*

baudrate: int = 9600

Baudrate for MBW973 is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, [hvl_ccb.comm.serial.SerialCommunicationParity](#)] = 'N'

MBW973 does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, [hvl_ccb.comm.serial.SerialCommunicationStopbits](#)] = 1

MBW973 does use one stop bit


terminator: bytes = b'\r'

The terminator is only CR

timeout: Union[int, float] = 3

use 3 seconds timeout as default

[hvl_ccb.dev.newport](#)

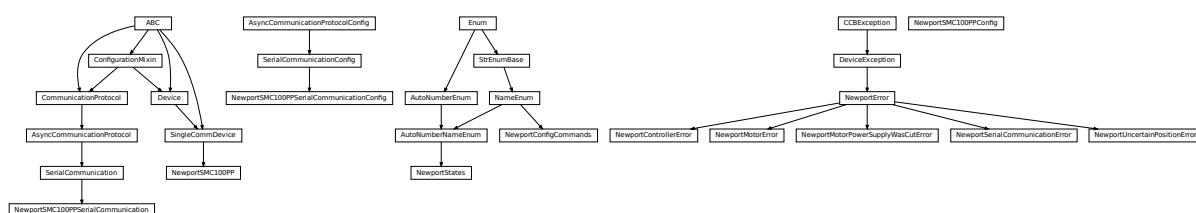

Device class for Newport SMC100PP stepper motor controller with serial communication.

The SMC100PP is a single axis motion controller/driver for stepper motors up to 48 VDC at 1.5 A rms. Up to 31 controllers can be networked through the internal RS-485 communication link.

Manufacturer homepage: <https://www.newport.com/f/sm100-single-axis-dc-or-stepper-motion-controller>

class NewportConfigCommands(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*, *boundary=None*)

Bases: [hvl_ccb.utils.enum.NameEnum](#)

Commands predefined by the communication protocol of the SMC100PP

AC = 'acceleration'

```

BA = 'backlash_compensation'
BH = 'hysteresis_compensation'
FRM = 'micro_step_per_full_step_factor'
FRS = 'motion_distance_per_full_step'
HT = 'home_search_type'
JR = 'jerk_time'
OH = 'home_search_velocity'
OT = 'home_search_timeout'
QIL = 'peak_output_current_limit'
SA = 'rs485_address'
SL = 'negative_software_limit'
SR = 'positive_software_limit'
VA = 'velocity'
VB = 'base_velocity'
ZX = 'stage_configuration'

exception NewportControllerError
    Bases: hvl\_ccb.dev.newport.NewportError
    Error with the Newport controller.

exception NewportError
    Bases: hvl\_ccb.dev.base.DeviceException
    General Exception for Newport Device

exception NewportMotorError
    Bases: hvl\_ccb.dev.newport.NewportError
    Error with the Newport motor.

exception NewportMotorPowerSupplyWasCutError
    Bases: hvl\_ccb.dev.newport.NewportError
    Error with the Newport motor after the power supply was cut and then restored, without interrupting the communication with the controller.

class NewportSMC100PP(com, dev_config=None)
    Bases: hvl\_ccb.dev.base.SingleCommDevice
    Device class of the Newport motor controller SMC100PP

class MotorErrors(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: aenum.Enum
    Possible motor errors reported by the motor during get_state().

    DC_VOLTAGE_TOO_LOW = 3
    FOLLOWING_ERROR = 6
    HOMING_TIMEOUT = 5
    NED_END_OF_TURN = 11

```

```
OUTPUT_POWER_EXCEEDED = 2
PEAK_CURRENT_LIMIT = 9
POS_END_OF_TURN = 10
RMS_CURRENT_LIMIT = 8
SHORT_CIRCUIT = 7
WRONG_ESP_STAGE = 4

class StateMessages(value=<no_arg>, names=None, module=None, type=None, start=1,
                     boundary=None)
Bases: aenum.Enum

Possible messages returned by the controller on get_state() query.

CONFIG = '14'
DISABLE_FROM_JOGGING = '3E'
DISABLE_FROM_MOVING = '3D'
DISABLE_FROM_READY = '3C'
HOMING_FROM_RS232 = '1E'
HOMING_FROM_SMC = '1F'
JOGGING_FROM_DISABLE = '47'
JOGGING_FROM_READY = '46'
MOVING = '28'
NO_REF_ESP_STAGE_ERROR = '10'
NO_REF_FROM_CONFIG = '0C'
NO_REF_FROM_DISABLED = '0D'
NO_REF_FROM_HOMING = '0B'
NO_REF_FROM_JOGGING = '11'
NO_REF_FROM_MOVING = '0F'
NO_REF_FROM_READY = '0E'
NO_REF_FROM_RESET = '0A'
READY_FROM_DISABLE = '34'
READY_FROM_HOMING = '32'
READY_FROM_JOGGING = '35'
READY_FROM_MOVING = '33'

States
alias of hvl\_ccb.dev.newport.NewportStates

static config_cls()
    Return the default configdataclass class.

        Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.
```

Returns the type of the standard communication protocol for this device

exit_configuration(*add*: *Optional[int]* = *None*) → *None*

Exit the CONFIGURATION state and go back to the NOT REFERENCED state. All configuration parameters are saved to the device's memory.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_acceleration(*add*: *Optional[int]* = *None*) → *Union[int, float]*

Leave the configuration state. The configuration parameters are saved to the device's memory.

Parameters **add** – controller address (1 to 31)

Returns acceleration (preset units/s^2), value between 1e-6 and 1e12

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_controller_information(*add*: *Optional[int]* = *None*) → *str*

Get information on the controller name and driver version

Parameters **add** – controller address (1 to 31)

Returns controller information

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_motor_configuration(*add*: *Optional[int]* = *None*) → *Dict[str, float]*

Query the motor configuration and returns it in a dictionary.

Parameters **add** – controller address (1 to 31)

Returns dictionary containing the motor's configuration

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

get_move_duration(*dist*: *Union[int, float]*, *add*: *Optional[int]* = *None*) → *float*

Estimate the time necessary to move the motor of the specified distance.

Parameters

- **dist** – distance to travel
- **add** – controller address (1 to 31), defaults to self.address

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_negative_software_limit(*add: Optional[int] = None*) → Union[int, float]

Get the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

Parameters *add* – controller address (1 to 31)

Returns negative software limit (preset units), value between -1e12 and 0

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_position(*add: Optional[int] = None*) → float

Returns the value of the current position.

Parameters *add* – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportUncertainPositionError* – if the position is ambiguous

get_positive_software_limit(*add: Optional[int] = None*) → Union[int, float]

Get the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

Parameters *add* – controller address (1 to 31)

Returns positive software limit (preset units), value between 0 and 1e12

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error

get_state(*add: int = None*) → *StateMessages*

Check on the motor errors and the controller state

Parameters *add* – controller address (1 to 31)

Raises

- *SerialCommunicationIOError* – if the com is closed
- *NewportSerialCommunicationError* – if an unexpected answer is obtained
- *NewportControllerError* – if the controller reports an error
- *NewportMotorError* – if the motor reports an error

Returns state message from the device (member of StateMessages)

go_home(*add: Optional[int] = None*) → None

Move the motor to its home position.

Parameters **add** – controller address (1 to 31), defaults to self.address

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

go_to_configuration(*add: Optional[int] = None*) → None

This method is executed during start(). It can also be executed after a reset(). The controller is put in CONFIG state, where configuration parameters can be changed.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

initialize(*add: Optional[int] = None*) → None

Puts the controller from the NOT_REF state to the READY state. Sends the motor to its “home” position.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

move_to_absolute_position(*pos: Union[int, float], add: Optional[int] = None*) → None

Move the motor to the specified position.

Parameters

- **pos** – target absolute position (affected by the configured offset)
- **add** – controller address (1 to 31), defaults to self.address

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

move_to_relative_position(*pos: Union[int, float], add: Optional[int] = None*) → None

Move the motor of the specified distance.

Parameters

- **pos** – distance to travel (the sign gives the direction)
- **add** – controller address (1 to 31), defaults to self.address

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

reset(*add*: *Optional[int]* = *None*) → *None*

Resets the controller, equivalent to a power-up. This puts the controller back to NOT REFERENCED state, which is necessary for configuring the controller.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_acceleration(*acc*: *Union[int, float]*, *add*: *Optional[int]* = *None*) → *None*

Leave the configuration state. The configuration parameters are saved to the device's memory.

Parameters

- **acc** – acceleration (preset units/s^2), value between 1e-6 and 1e12
- **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_motor_configuration(*add*: *Optional[int]* = *None*, *config*: *Optional[dict]* = *None*) → *None*

Set the motor configuration. The motor must be in CONFIG state.

Parameters

- **add** – controller address (1 to 31)
- **config** – dictionary containing the motor's configuration

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

set_negative_software_limit(*lim*: *Union[int, float]*, *add*: *Optional[int]* = *None*) → *None*

Set the negative software limit (the maximum position that the motor is allowed to travel to towards the left).

Parameters

- **lim** – negative software limit (preset units), value between -1e12 and 0
- **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained

- **NewportControllerError** – if the controller reports an error

set_positive_software_limit(*lim: Union[int, float], add: Optional[int] = None*) → None

Set the positive software limit (the maximum position that the motor is allowed to travel to towards the right).

Parameters

- **lim** – positive software limit (preset units), value between 0 and 1e12
- **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

start()

Opens the communication protocol and applies the config.

Raises **SerialCommunicationIOError** – when communication port cannot be opened

stop() → None

Stop the device. Close the communication protocol.

stop_motion(*add: Optional[int] = None*) → None

Stop a move in progress by decelerating the positioner immediately with the configured acceleration until it stops. If a controller address is provided, stops a move in progress on this controller, else stops the moves on all controllers.

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

wait_until_motor_initialized(*add: Optional[int] = None*) → None

Wait until the motor leaves the HOMING state (at which point it should have arrived to the home position).

Parameters **add** – controller address (1 to 31)

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

```
class NewportSMC100PPConfig(address: int = 1, user_position_offset: Union[int, float] = 23.987,
    screw_scaling: Union[int, float] = 1, exit_configuration_wait_sec: Union[int,
    float] = 5, move_wait_sec: Union[int, float] = 1, acceleration: Union[int, float]
    = 10, backlash_compensation: Union[int, float] = 0, hysteresis_compensation:
    Union[int, float] = 0.015, micro_step_per_full_step_factor: int = 100,
    motion_distance_per_full_step: Union[int, float] = 0.01, home_search_type:
    Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] =
    HomeSearch.HomeSwitch, jerk_time: Union[int, float] = 0.04,
    home_search_velocity: Union[int, float] = 4, home_search_timeout: Union[int,
    float] = 27.5, home_search_polling_interval: Union[int, float] = 1,
    peak_output_current_limit: Union[int, float] = 0.4, rs485_address: int = 2,
    negative_software_limit: Union[int, float] = - 23.5, positive_software_limit:
    Union[int, float] = 25, velocity: Union[int, float] = 4, base_velocity: Union[int,
    float] = 0, stage_configuration: Union[int,
    hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageConfig] =
    EspStageConfig.EnableEspStageCheck)
```

Bases: `object`

Configuration dataclass for the Newport motor controller SMC100PP.

```
class EspStageConfig(value=<no_arg>, names=None, module=None, type=None, start=1,
    boundary=None)
```

Bases: `aenum.IntEnum`

Different configurations to check or not the motor configuration upon power-up.

`DisableEspStageCheck` = 1

`EnableEspStageCheck` = 3

`UpdateEspStageInfo` = 2

```
class HomeSearch(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
```

Bases: `aenum.IntEnum`

Different methods for the motor to search its home position during initialization.

`CurrentPosition` = 1

`EndOfRunSwitch` = 4

`EndOfRunSwitch_and_Index` = 3

`HomeSwitch` = 2

`HomeSwitch_and_Index` = 0

`acceleration: Union[int, float]` = 10

`address: int` = 1

`backlash_compensation: Union[int, float]` = 0

`base_velocity: Union[int, float]` = 0

`clean_values()`

`exit_configuration_wait_sec: Union[int, float]` = 5

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

home_search_polling_interval: Union[int, float] = 1

home_search_timeout: Union[int, float] = 27.5

home_search_type: Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.HomeSearch] = 2

home_search_velocity: Union[int, float] = 4

hysteresis_compensation: Union[int, float] = 0.015

is_configdataclass = True

jerk_time: Union[int, float] = 0.04

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

micro_step_per_full_step_factor: int = 100

motion_distance_per_full_step: Union[int, float] = 0.01

property motor_config: Dict[str, float]

Gather the configuration parameters of the motor into a dictionary.

Returns dict containing the configuration parameters of the motor

move_wait_sec: Union[int, float] = 1

negative_software_limit: Union[int, float] = -23.5

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

peak_output_current_limit: Union[int, float] = 0.4

positive_software_limit: Union[int, float] = 25

post_force_value(fieldname, value)

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

rs485_address: int = 2

screw_scaling: Union[int, float] = 1

stage_configuration: Union[int, hvl_ccb.dev.newport.NewportSMC100PPConfig.EspStageConfig] = 3

user_position_offset: Union[int, float] = 23.987

velocity: Union[int, float] = 4

```
class NewportSMC100PPSerialCommunication(configuration)
```

Bases: [hvl_ccb.comm.serial.SerialCommunication](#)

Specific communication protocol implementation for NewportSMC100 controller. Already predefines device-specific protocol parameters in config.

```
class ControllerErrors(value=<no_arg>, names=None, module=None, type=None, start=1,  
boundary=None)
```

Bases: [aenum.Enum](#)

Possible controller errors with values as returned by the device in response to sent commands.

ADDR_INCORRECT = 'B'

CMD_EXEC_ERROR = 'V'

CMD_NOT_ALLOWED = 'D'

CMD_NOT_ALLOWED_CC = 'X'

CMD_NOT_ALLOWED_CONFIGURATION = 'I'

CMD_NOT_ALLOWED_DISABLE = 'J'

CMD_NOT_ALLOWED_HOMING = 'L'

CMD_NOT_ALLOWED_MOVING = 'M'

CMD_NOT_ALLOWED_NOT_REFERENCED = 'H'

CMD_NOT_ALLOWED_PP = 'W'

CMD_NOT_ALLOWED_READY = 'K'

CODE_OR_ADDR_INVALID = 'A'

COM_TIMEOUT = 'S'

DISPLACEMENT_OUT_OF_LIMIT = 'G'

EEPROM_ACCESS_ERROR = 'U'

ESP_STAGE_NAME_INVALID = 'F'

HOME_STARTED = 'E'

NO_ERROR = '@'

PARAM_MISSING_OR_INVALID = 'C'

POSITION_OUT_OF_LIMIT = 'N'

check_for_error(add: int) → None

Ask the Newport controller for the last error it recorded.

This method is called after every command or query.

Parameters add – controller address (1 to 31)

Raises

- [SerialCommunicationIOError](#) – if the com is closed
- [NewportSerialCommunicationError](#) – if an unexpected answer is obtained
- [NewportControllerError](#) – if the controller reports an error

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(*add*: int, *cmd*: str, *param*: Optional[Union[int, float, str]] = None) → str

Send a query to the controller, read the answer, and check for errors. The prefix add+cmd is removed from the answer.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **param** – optional parameter (int/float/str) appended to the command

Returns the answer from the device without the prefix

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

query_multiple(*add*: int, *cmd*: str, *prefixes*: List[str]) → List[str]

Send a query to the controller, read the answers, and check for errors. The prefixes are removed from the answers.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent
- **prefixes** – prefixes of each line expected in the answer

Returns list of answers from the device without prefix

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

read_text() → str

Read one line of text from the serial port, and check for presence of a null char which indicates that the motor power supply was cut and then restored. The input buffer may hold additional data afterwards, since only one line is read.

This method uses *self.access_lock* to ensure thread-safety.

Returns String read from the serial port; '' if there was nothing to read.

Raises

- **SerialCommunicationIOError** – when communication port is not opened
- **NewportMotorPowerSupplyWasCutError** – if a null char is read

send_command(*add*: int, *cmd*: str, *param*: Optional[Union[int, float, str]] = None) → None

Send a command to the controller, and check for errors.

Parameters

- **add** – the controller address (1 to 31)
- **cmd** – the command to be sent

- **param** – optional parameter (int/float/str) appended to the command

Raises

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained
- **NewportControllerError** – if the controller reports an error

send_stop(*add*: int) → None

Send the general stop ST command to the controller, and check for errors.

Parameters *add* – the controller address (1 to 31)**Returns** ControllerErrors reported by Newport Controller**Raises**

- **SerialCommunicationIOError** – if the com is closed
- **NewportSerialCommunicationError** – if an unexpected answer is obtained

```
class NewportSMC100PPSSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'ascii',
                                                encoding_error_handling: str = 'replace',
                                                wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                                default_n_attempts_read_text_nonempty: int = 10,
                                                port: Union[str, NoneType] = None, baudrate: int =
                                                57600, parity: Union[str,
                                                hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                Union[int,
                                                hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                <SerialCommunicationBytesize.EIGHTBITS: 8>,
                                                timeout: Union[int, float] = 10)
```

Bases: *hvl_ccb.comm.serial.SerialCommunicationConfig***baudrate: int = 57600**

Baudrate for NewportSMC100 controller is 57600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

NewportSMC100 controller uses 8 bits for one data byte

encoding: str = 'ascii'

use ASCII as de-/encoding, cf. the manual

encoding_error_handling: str = 'replace'

replace bytes with instead of raising utf-8 exception when decoding fails

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.**Parameters**

- **fieldname** – name of the field
- **value** – value to assign

```
classmethod keys() → Sequence[str]
    Returns a list of all configdataclass fields key-names.

    Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]
    Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

    Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'
    NewportSMC100 controller does not use parity

classmethod required_keys() → Sequence[str]
    Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

    Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
    NewportSMC100 controller uses one stop bit

terminator: bytes = b'\r\n'
    The terminator is CR/LF

timeout: Union[int, float] = 10
    use 10 seconds timeout as default

exception NewportSerialCommunicationError
    Bases: hvl_ccb.dev.newport.NewportError
    Communication error with the Newport controller.


class NewportStates(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)
    Bases: hvl_ccb.utils.enum.AutoNumberNameEnum

    States of the Newport controller. Certain commands are allowed only in certain states.

    CONFIG = 3
    DISABLE = 6
    HOMING = 2
    JOGGING = 7
    MOVING = 5
    NO_REF = 1
    READY = 4

exception NewportUncertainPositionError
    Bases: hvl_ccb.dev.newport.NewportError
    Error with the position of the Newport motor.
```

`hvl_ccb.dev.pfeiffer_tpg`

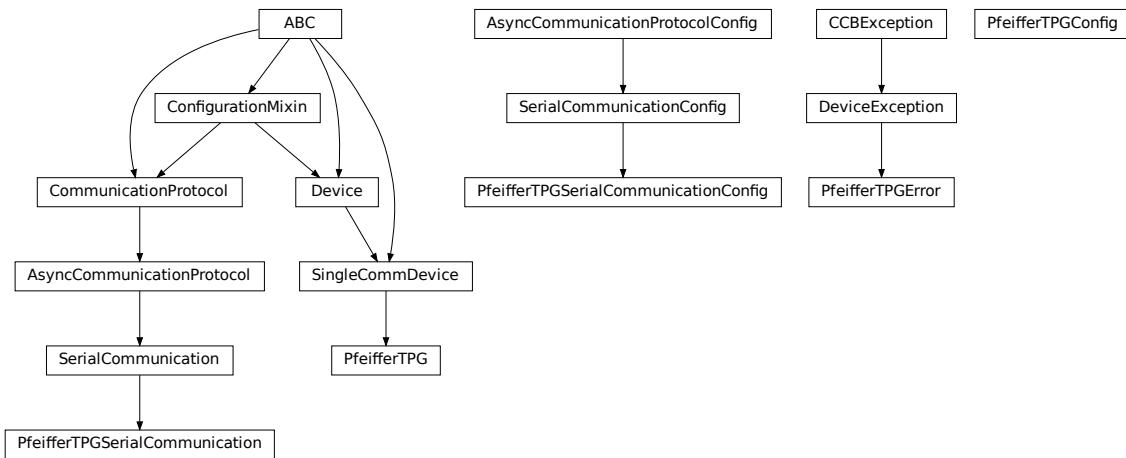

Device class for Pfeiffer TPG controllers.

The Pfeiffer TPG control units are used to control Pfeiffer Compact Gauges. Models: TPG 251 A, TPG 252 A, TPG 256A, TPG 261, TPG 262, TPG 361, TPG 362 and TPG 366.

Manufacturer homepage: <https://www.pfeiffer-vacuum.com/en/products/measurement-analysis/> measurement/activeline/controllers/

class PfeifferTPG(*com, dev_config=None*)
Bases: *hvl_ccb.dev.base.SingleCommDevice*

Pfeiffer TPG control unit device class

class PressureUnits(*value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None*)
Bases: *hvl_ccb.utils.enum.NameEnum*

Enum of available pressure units for the digital display. “0” corresponds either to bar or to mbar depending on the TPG model. In case of doubt, the unit is visible on the digital display.

Micron = 3

Pascal = 2

Torr = 1

Volt = 5

bar = 0

hPascal = 4

mbar = 0

class SensorStatus(*value*)

Bases: *enum.IntEnum*

An enumeration.

Identification_error = 6

No_sensor = 5

```

Ok = 0
OVERRANGE = 2
Sensor_error = 3
Sensor_off = 4
Underrange = 1

class SensorTypes(value)
    Bases: enum.Enum

    An enumeration.

    CMR = 4
    IKR = 2
    IKR11 = 2
    IKR9 = 2
    IMR = 5
    None = 7
    PBR = 6
    PKR = 3
    TPR = 1
    noSENSOR = 7
    noSen = 7

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

get_full_scale_mbar() → List[Union[int, float]]
    Get the full scale range of the attached sensors

    Returns full scale range values in mbar, like [0.01, 1, 0.1, 1000, 50000, 10]

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • PfeifferTPGError – if command fails

get_full_scale_unitless() → List[int]
    Get the full scale range of the attached sensors. See lookup table between command and corresponding
    pressure in the device user manual.

    Returns list of full scale range values, like [0, 1, 3, 3, 2, 0]

    Raises
        • SerialCommunicationIOError – when communication port is not opened
        • PfeifferTPGError – if command fails

```

identify_sensors() → None

Send identification request TID to sensors on all channels.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

measure(channel: int) → Tuple[str, float]

Get the status and measurement of one sensor

Parameters **channel** – int channel on which the sensor is connected, with $1 \leq \text{channel} \leq \text{number_of_sensors}$

Returns measured value as float if measurement successful, sensor status as string if not

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

measure_all() → List[Tuple[str, float]]

Get the status and measurement of all sensors (this command is not available on all models)

Returns list of measured values as float if measurements successful, and or sensor status as strings if not

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

property number_of_sensors

set_display_unit(unit: Union[str, hvl_ccb.dev.pfeiffer_tpg.PfeifferTPG.PressureUnits]) → None

Set the unit in which the measurements are shown on the display.

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

set_full_scale_mbar(fsr: List[Union[int, float]]) → None

Set the full scale range of the attached sensors (in unit mbar)

Parameters **fsr** – full scale range values in mbar, for example [0.01, 1000]

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

set_full_scale_unitless(fsr: List[int]) → None

Set the full scale range of the attached sensors. See lookup table between command and corresponding pressure in the device user manual.

Parameters **fsr** – list of full scale range values, like [0, 1, 3, 3, 2, 0]

Raises

- *SerialCommunicationIOError* – when communication port is not opened
- *PfeifferTPGError* – if command fails

start() → None

Start this device. Opens the communication protocol, and identify the sensors.

Raises `SerialCommunicationIOError` – when communication port cannot be opened

stop() → None

Stop the device. Closes also the communication protocol.

property unit

The pressure unit of readings is always mbar, regardless of the display unit.

class PfeifferTPGConfig(*model*: Union[str, `hvl_ccb.dev.pfeiffer_tpg.PfeifferTPGConfig.Model`] = `Model.TPG25xA`)

Bases: object

Device configuration dataclass for Pfeiffer TPG controllers.

class Model(*value*=<no_arg>, *names*=None, *module*=None, *type*=None, *start*=1, *boundary*=None)

Bases: `hvl_ccb.utils.enum.NameEnum`

An enumeration.

`TPG25xA` = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

`TPGx6x` = {0.01: 0, 0.1: 1, 1: 2, 10: 3, 100: 4, 1000: 5, 2000: 6, 5000: 7, 10000: 8, 50000: 9}

is_valid_scale_range_reversed_str(*v*: str) → bool

Check if given string represents a valid reversed scale range of a model.

Parameters *v* – Reversed scale range string.

Returns *True* if valid, *False* otherwise.

clean_values()

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

model: Union[str, `hvl_ccb.dev.pfeiffer_tpg.PfeifferTPGConfig.Model`] = {0.1: 8, 1: 0, 10: 1, 100: 2, 1000: 3, 2000: 4, 5000: 5, 10000: 6, 50000: 7}

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception PfeifferTPGError

Bases: *hvl_ccb.dev.base.DeviceException*

Error with the Pfeiffer TPG Controller.

class PfeifferTPGSerialCommunication(configuration)

Bases: *hvl_ccb.comm.serial.SerialCommunication*

Specific communication protocol implementation for Pfeiffer TPG controllers. Already predefines device-specific protocol parameters in config.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

query(cmd: str) → str

Send a query, then read and returns the first line from the com port.

Parameters cmd – query message to send to the device

Returns first line read on the com

Raises

- **SerialCommunicationIOError** – when communication port is not opened

- **PfeifferTPGError** – if the device does not acknowledge the command or if the answer from the device is empty

send_command(cmd: str) → None

Send a command to the device and check for acknowledgement.

Parameters cmd – command to send to the device

Raises

- **SerialCommunicationIOError** – when communication port is not opened

- **PfeifferTPGError** – if the answer from the device differs from the expected acknowledgement character ‘chr(6)’.

```
class PfeifferTPGSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8',
                                             encoding_error_handling: str = 'strict',
                                             wait_sec_read_text_nonempty: Union[int, float] = 0.5,
                                             default_n_attempts_read_text_nonempty: int = 10, port:
                                             Union[str, NoneType] = None, baudrate: int = 9600, parity:
                                             Union[str,
                                                   hvl_ccb.comm.serial.SerialCommunicationParity] =
                                                   <SerialCommunicationParity.NONE: 'N'>, stopbits:
                                                   Union[int,
                                                       hvl_ccb.comm.serial.SerialCommunicationStopbits] =
                                                       <SerialCommunicationStopbits.ONE: 1>, bytesize:
                                                       Union[int,
                                                       hvl_ccb.comm.serial.SerialCommunicationBytesize] =
                                                       <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout:
                                                       Union[int, float] = 3)
```

Bases: `hvl_ccb.comm.serial.SerialCommunicationConfig`

baudrate: int = 9600

Baudrate for Pfeiffer TPG controllers is 9600 baud

bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8

One byte is eight bits long

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'

Pfeiffer TPG controllers do not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1

Pfeiffer TPG controllers use one stop bit

terminator: bytes = b'\r\n'

The terminator is <CR><LF>

```
timeout: Union[int, float] = 3
use 3 seconds timeout as default
```

`hvl_ccb.dev.picotech_pt104`


NOTE: PicoSDK Python wrappers already on import attempt to load the PicoSDK library; thus, the API docs can only be generated in a system with the latter installed and are by default disabled.

To build the API docs for this submodule locally edit the `docs/hvl_ccb.dev.picotech_pt104.rst` file to remove the `.. code-block::` directive preceding the following directives:

```
.. inheritance-diagram:: hvl_ccb.dev.picotech_pt104
:parts: 1

.. automodule:: hvl_ccb.dev.picotech_pt104
:members:
:undoc-members:
:show-inheritance:
```

`hvl_ccb.dev.rs_rto1024`

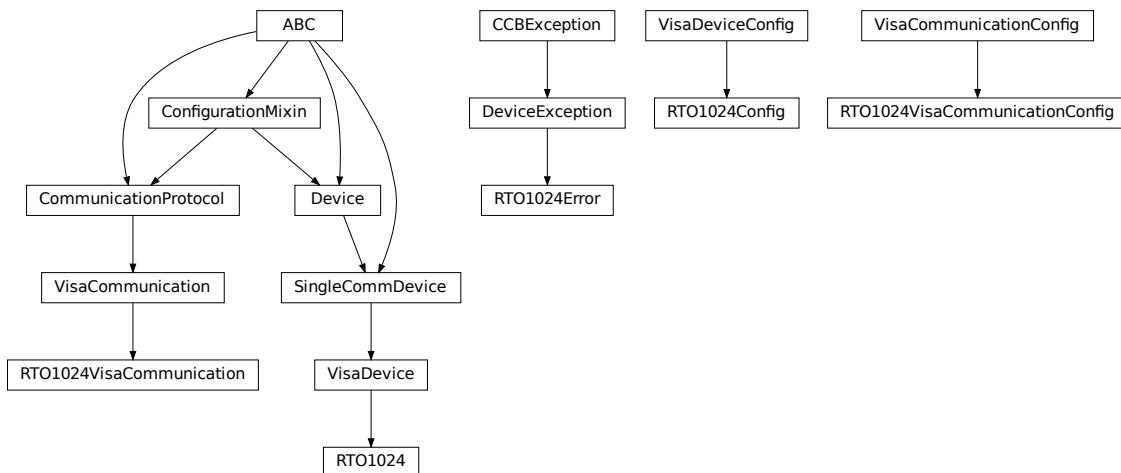

Python module for the Rhode & Schwarz RTO 1024 oscilloscope. The communication to the device is through VISA, type TCPIP / INSTR.

```
class RT01024(com: Union[hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunication,
                        hvl_ccb.dev.rs_rto1024.RTO1024VisaCommunicationConfig, dict], dev_config:
                        Union[hvl_ccb.dev.rs_rto1024.RTO1024Config, dict])
Bases: hvl_ccb.dev.visa.VisaDevice
```

Device class for the Rhode & Schwarz RTO 1024 oscilloscope.

```
class TriggerModes(value=<no_arg>, names=None, module=None, type=None, start=1,
                   boundary=None)
Bases: hvl_ccb.utils.enum.AutoNumberNameEnum
```

Enumeration for the three available trigger modes.

AUTO = 1

FREERUN = 3

NORMAL = 2

classmethod names()

Returns a list of the available trigger modes. :return: list of strings

activate_measurements(meas_n: int, source: str, measurements: List[str], category: str = 'AMPTime')

Activate the list of ‘measurements’ of the waveform ‘source’ in the measurement box number ‘meas_n’. The list ‘measurements’ starts with the main measurement and continues with additional measurements of the same ‘category’.

Parameters

- **meas_n** – measurement number 1..8
- **source** – measurement source, for example C1W1
- **measurements** – list of measurements, the first one will be the main measurement.
- **category** – the category of measurements, by default AMPTime

backup_waveform(filename: str) → None

Backup a waveform file from the standard directory specified in the device configuration to the standard backup destination specified in the device configuration. The filename has to be specified without .bin or path.

Parameters filename – The waveform filename without extension and path

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Return the default communication protocol for this device type, which is VisaCommunication.

Returns the VisaCommunication class

file_copy(source: str, destination: str) → None

Copy a file from one destination to another on the oscilloscope drive. If the destination file already exists, it is overwritten without notice.

Parameters

- **source** – absolute path to the source file on the DSO filesystem
- **destination** – absolute path to the destination file on the DSO filesystem

Raises RT01024Error – if the operation did not complete

get_acquire_length() → float

Gets the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]

Returns the time for one acquisition. Range: 250e-12 ... 500 [s]

get_channel_offset(channel: int) → float

Gets the voltage offset of the indicated channel.

Parameters `channel` – is the channel number (1..4)

Returns channel offset voltage in V (value between -1 and 1)

get_channel_position(`channel: int`) → float

Gets the vertical position of the indicated channel.

Parameters `channel` – is the channel number (1..4)

Returns channel position in div (value between -5 and 5)

get_channel_range(`channel: int`) → float

Queries the channel range in V.

Parameters `channel` – is the input channel (1..4)

Returns channel range in V

get_channel_scale(`channel: int`) → float

Queries the channel scale in V/div.

Parameters `channel` – is the input channel (1..4)

Returns channel scale in V/div

get_channel_state(`channel: int`) → bool

Queries if the channel is active or not.

Parameters `channel` – is the input channel (1..4)

Returns True if active, else False

get_reference_point() → int

Gets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]

Returns the reference in %

get_repetitions() → int

Get the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- *RST = 1

Returns the number of waveforms to acquire

get_timestamps() → List[float]

Gets the timestamps of all recorded frames in the history and returns them as a list of floats.

Returns list of timestamps in [s]

Raises `RT01024Error` – if the timestamps are invalid

list_directory(`path: str`) → List[Tuple[str, str, int]]

List the contents of a given directory on the oscilloscope filesystem.

Parameters `path` – is the path to a folder

Returns a list of filenames in the given folder

load_configuration(*filename*: str) → None

Load current settings from a configuration file. The filename has to be specified without base directory and ‘.dfl’ extension.

Information from the manual *ReCaLL* calls up the instrument settings from an intermediate memory identified by the specified number. The instrument settings can be stored to this memory using the command *SAV with the associated number. It also activates the instrument settings which are stored in a file and loaded using *MMEMemory:LOAD:STATE*.

Parameters **filename** – is the name of the settings file without path and extension

local_display(*state*: bool) → None

Enable or disable local display of the scope.

Parameters **state** – is the desired local display state

prepare_ultra_segmentation() → None

Make ready for a new acquisition in ultra segmentation mode. This function does one acquisition without ultra segmentation to clear the history and prepare for a new measurement.

read_measurement(*meas_n*: int, *name*: str) → float

Parameters

- **meas_n** – measurement number 1..8
- **name** – measurement name, for example “MAX”

Returns measured value

run_continuous_acquisition() → None

Start acquiring continuously.

run_single_acquisition() → None

Start a single or Nx acquisition.

save_configuration(*filename*: str) → None

Save the current oscilloscope settings to a file. The filename has to be specified without path and ‘.dfl’ extension, the file will be saved to the configured settings directory.

Information from the manual *SAVe* stores the current instrument settings under the specified number in an intermediate memory. The settings can be recalled using the command *RCL with the associated number. To transfer the stored instrument settings to a file, use *MMEMemory:STORe:STATE*.

Parameters **filename** – is the name of the settings file without path and extension

save_waveform_history(*filename*: str, *channel*: int, *waveform*: int = 1) → None

Save the history of one channel and one waveform to a .bin file. This function is used after an acquisition using sequence trigger mode (with or without ultra segmentation) was performed.

Parameters

- **filename** – is the name (without extension) of the file
- **channel** – is the channel number
- **waveform** – is the waveform number (typically 1)

Raises *RT01024Error* – if storing waveform times out

set_acquire_length(*timerange*: float) → None

Defines the time of one acquisition, that is the time across the 10 divisions of the diagram.

- Range: 250E-12 ... 500 [s]
- Increment: 1E-12 [s]
- *RST = 0.5 [s]

Parameters **timerange** – is the time for one acquisition. Range: 250e-12 ... 500 [s]

set_channel_offset(*channel: int, offset: float*) → None

Sets the voltage offset of the indicated channel.

- Range: Dependent on the channel scale and coupling [V]
- Increment: Minimum 0.001 [V], may be higher depending on the channel scale and coupling
- *RST = 0

Parameters

- **channel** – is the channel number (1..4)
- **offset** – Offset voltage. Positive values move the waveform down, negative values move it up.

set_channel_position(*channel: int, position: float*) → None

Sets the vertical position of the indicated channel as a graphical value.

- Range: -5.0 ... 5.0 [div]
- Increment: 0.02
- *RST = 0

Parameters

- **channel** – is the channel number (1..4)
- **position** – is the position. Positive values move the waveform up, negative values move it down.

set_channel_range(*channel: int, v_range: float*) → None

Sets the voltage range across the 10 vertical divisions of the diagram. Use the command alternatively instead of set_channel_scale.

- Range for range: Depends on attenuation factors and coupling. With 1:1 probe and external attenuations and 50 input coupling, the range is 10 mV to 10 V. For 1 M input coupling, it is 10 mV to 100 V. If the probe and/or external attenuation is changed, multiply the range values by the attenuation factors.
- Increment: 0.01
- *RST = 0.5

Parameters

- **channel** – is the channel number (1..4)
- **v_range** – is the vertical range [V]

set_channel_scale(*channel: int, scale: float*) → None

Sets the vertical scale for the indicated channel. The scale value is given in volts per division.

- Range for scale: depends on attenuation factor and coupling. With 1:1 probe and external attenuations and 50 input coupling, the vertical scale (input sensitivity) is 1 mV/div to 1 V/div. For 1 M input coupling, it is 1 mV/div to 10 V/div. If the probe and/or external attenuation is changed, multiply the values by the attenuation factors to get the actual scale range.
- Increment: 1e-3
- *RST = 0.05

See also: `set_channel_range`

Parameters

- **channel** – is the channel number (1..4)
- **scale** – is the vertical scaling [V/div]

set_channel_state(*channel: int, state: bool*) → None

Switches the channel signal on or off.

Parameters

- **channel** – is the input channel (1..4)
- **state** – is True for on, False for off

set_reference_point(*percentage: int*) → None

Sets the reference point of the time scale in % of the display. If the “Trigger offset” is zero, the trigger point matches the reference point. ReferencePoint = zero pint of the time scale

- Range: 0 … 100 [%]
- Increment: 1 [%]
- *RST = 50 [%]

Parameters **percentage** – is the reference in %

set_repetitions(*number: int*) → None

Set the number of acquired waveforms with RUN Nx SINGLE. Also defines the number of waveforms used to calculate the average waveform.

- Range: 1 … 16777215
- Increment: 10
- *RST = 1

Parameters **number** – is the number of waveforms to acquire

set_trigger_level(*channel: int, level: float, event_type: int = 1*) → None

Sets the trigger level for the specified event and source.

- Range: -10 to 10 V
- Increment: 1e-3 V
- *RST = 0 V

Parameters

- **channel** – indicates the trigger source.
 - 1..4 = channel 1 to 4, available for all event types 1..3

- 5 = external trigger input on the rear panel for analog signals, available for A-event type
= 1
- 6..9 = not available
- **level** – is the voltage for the trigger level in [V].
- **event_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

set_trigger_mode(*mode*: Union[str, hvl_ccb.dev.rs_rto1024.RTO1024.TriggerModes]) → None
Sets the trigger mode which determines the behavior of the instrument if no trigger occurs.

Parameters **mode** – is either auto, normal, or freerun.

Raises **RT01024Error** – if an invalid triggermode is selected

set_trigger_source(*channel*: int, *event_type*: int = 1) → None
Set the trigger (Event A) source channel.

Parameters

- **channel** – is the channel number (1..4)
- **event_type** – is the event type. 1: A-Event, 2: B-Event, 3: R-Event

start() → None

Start the RTO1024 oscilloscope and bring it into a defined state and remote mode.

stop() → None

Stop the RTO1024 oscilloscope, reset events and close communication. Brings back the device to a state where local operation is possible.

stop_acquisition() → None
Stop any acquisition.

class RT01024Config(*waveforms_path*: str, *settings_path*: str, *backup_path*: str, *spoll_interval*: Union[int, float] = 0.5, *spoll_start_delay*: Union[int, float] = 2, *command_timeout_seconds*: Union[int, float] = 60, *wait_sec_short_pause*: Union[int, float] = 0.1, *wait_sec_enable_history*: Union[int, float] = 1, *wait_sec_post_acquisition_start*: Union[int, float] = 2)
Bases: [hvl_ccb.dev.visa.VisaDeviceConfig](#), [hvl_ccb.dev.rs_rto1024._RT01024ConfigDefaultsBase](#), [hvl_ccb.dev.rs_rto1024._RT01024ConfigBase](#)

Configdataclass for the RTO1024 device.

force_value(*fieldname*, *value*)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

exception RT01024Error

Bases: `hvl_ccb.dev.base.DeviceException`

class RT01024VisaCommunication(configuration)

Bases: `hvl_ccb.comm.visa.VisaCommunication`

Specialization of VisaCommunication for the RTO1024 oscilloscope

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class RT01024VisaCommunicationConfig(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], interface_type: Union[str, `hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType`] = InterfaceType.TCPIP_INSTR, board: int = 0, port: int = 5025, timeout: int = 5000, chunk_size: int = 204800, open_timeout: int = 1000, write_termination: str = '\n', read_termination: str = '\n', visa_backend: str = '')

Bases: `hvl_ccb.comm.visa.VisaCommunicationConfig`

Configuration dataclass for VisaCommunication with specifications for the RTO1024 device class.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after `value` has been forced on `fieldname`.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

interface_type: Union[str, `hvl_ccb.comm.visa.VisaCommunicationConfig.InterfaceType`] = 2

Interface type of the VISA connection, being one of `InterfaceType`.

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.


Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

[hvl_ccb.dev.se_il2t](#)

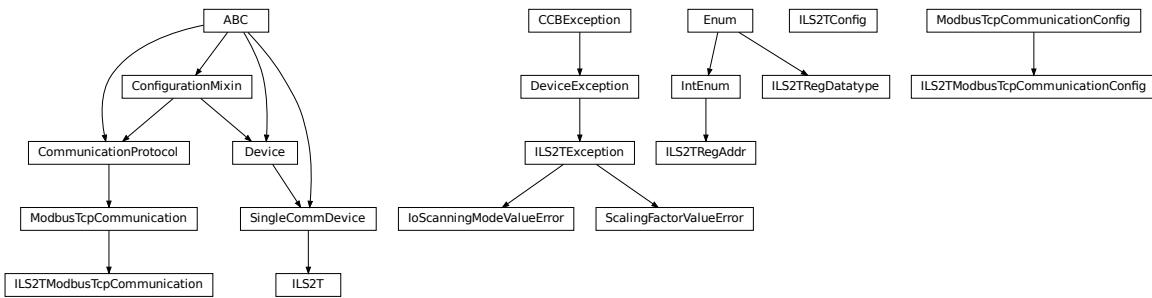

Device class for controlling a Schneider Electric ILS2T stepper drive over modbus TCP.

class ILS2T(com, dev_config=None)
Bases: [hvl_ccb.dev.base.SingleCommDevice](#)

Schneider Electric ILS2T stepper drive class.

ACTION_JOG_VALUE = 0

The single action value for *ILS2T.Mode.JOG*

class ActionsPtp(value)
Bases: [enum.IntEnum](#)

Allowed actions in the point to point mode (*ILS2T.Mode.PTP*).

ABSOLUTE_POSITION = 0

RELATIVE_POSITION_MOTOR = 2

RELATIVE_POSITION_TARGET = 1

DEFAULT_IO_SCANNING_CONTROL_VALUES = {'action': 2, 'continue_after_stop_cu': 0, 'disable_driver_di': 0, 'enable_driver_en': 0, 'execute_stop_sh': 0, 'fault_reset_fr': 0, 'mode': 3, 'quick_stop_qs': 0, 'ref_16': 1500, 'ref_32': 0, 'reset_stop_ch': 0}

Default IO Scanning control mode values

class Mode(value)
Bases: [enum.IntEnum](#)

ILS2T device modes

JOG = 1

PTP = 3

class Ref16Jog(value)
Bases: [enum.Flag](#)

Allowed values for ILS2T ref_16 register (the shown values are the integer representation of the bits), all in Jog mode = 1

FAST = 4

NEG = 2

NEG_FAST = 6

NONE = 0

```

POS = 1
POS_FAST = 5

RegAddr
    Modbus Register Adresses
    alias of hvl\_ccb.dev.se\_ilst2t.ILS2TRegAddr

RegDatatype
    Modbus Register Datatypes
    alias of hvl\_ccb.dev.se\_ilst2t.ILS2TRegDatatype

class State(value)
    Bases: enum.IntEnum
    State machine status values
    ON = 6
    QUICKSTOP = 7
    READY = 4

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls()
    Get the class for the default communication protocol used with this device.

    Returns the type of the standard communication protocol for this device

disable(log_warn: bool = True, wait_sec_max: Optional[int] = None) → bool
    Disable the driver of the stepper motor and enable the brake.

    Note: the driver cannot be disabled if the motor is still running.

Parameters

- log_warn – if log a warning in case the motor cannot be disabled.
- wait_sec_max – maximal wait time for the motor to stop running and to disable it; by default, with None, use a config value

Returns True if disable request could and was sent, False otherwise.

do_ioscanning_write(**kwargs: int) → None
    Perform a write operation using IO Scanning mode.

    Parameters kwargs – Keyword-argument list with options to send, remaining are taken from the defaults.

enable() → None
    Enable the driver of the stepper motor and disable the brake.

execute_absolute_position(position: int) → bool
    Execute a absolute position change, i.e. enable motor, perform absolute position change, wait until done and disable motor afterwards.

    Check position at the end if wrong do not raise error; instead just log and return check result.

    Parameters position – absolute position of motor in user defined steps.

    Returns True if actual position is as expected, False otherwise.

```

execute_relative_step(*steps: int*) → bool

Execute a relative step, i.e. enable motor, perform relative steps, wait until done and disable motor afterwards.

Check position at the end if wrong do not raise error; instead just log and return check result.

Parameters **steps** – Number of steps.

Returns *True* if actual position is as expected, *False* otherwise.

get_dc_volt() → float

Read the DC supply voltage of the motor.

Returns DC input voltage.

get_error_code() → Dict[int, Dict[str, Any]]

Read all messages in fault memory. Will read the full error message and return the decoded values. At the end the fault memory of the motor will be deleted. In addition, *reset_error* is called to re-enable the motor for operation.

Returns Dictionary with all information

get_position() → int

Read the position of the drive and store into status.

Returns Position step value

get_status() → Dict[str, int]

Perform an IO Scanning read and return the status of the motor.

Returns dict with status information.

get_temperature() → int

Read the temperature of the motor.

Returns Temperature in degrees Celsius.

jog_run(*direction: bool = True, fast: bool = False*) → None

Slowly turn the motor in positive direction.

jog_stop() → None

Stop turning the motor in Jog mode.

quickstop() → None

Stops the motor with high deceleration rate and falls into error state. Reset with *reset_error* to recover into normal state.

reset_error() → None

Resets the motor into normal state after quick stop or another error occurred.

set_jog_speed(*slow: int = 60, fast: int = 180*) → None

Set the speed for jog mode. Default values correspond to startup values of the motor.

Parameters

- **slow** – RPM for slow jog mode.
- **fast** – RPM for fast jog mode.

set_max_acceleration(*rpm_minute: int*) → None

Set the maximum acceleration of the motor.

Parameters **rpm_minute** – revolution per minute per minute

set_max_deceleration(*rpm_minute: int*) → None

Set the maximum deceleration of the motor.

Parameters `rpm_minute` – revolution per minute per minute

`set_max_rpm(rpm: int) → None`

Set the maximum RPM.

Parameters `rpm` – revolution per minute ($0 < rpm \leq RPM_MAX$)

Raises `ILS2TException` – if RPM is out of range

`set_ramp_type(ramp_type: int = -1) → None`

Set the ramp type. There are two options available: 0: linear ramp -1: motor optimized ramp

Parameters `ramp_type` – 0: linear ramp | -1: motor optimized ramp

`start() → None`

Start this device.

`stop() → None`

Stop this device. Disables the motor (applies brake), disables access and closes the communication protocol.

`user_steps(steps: int = 16384, revolutions: int = 1) → None`

Define steps per revolution. Default is 16384 steps per revolution. Maximum precision is 32768 steps per revolution.

Parameters

- **steps** – number of steps in *revolutions*.
- **revolutions** – number of revolutions corresponding to *steps*.

`write_absolute_position(position: int) → None`

Write instruction to turn the motor until it reaches the absolute position. This function does not enable or disable the motor automatically.

Parameters `position` – absolute position of motor in user defined steps.

`write_relative_step(steps: int) → None`

Write instruction to turn the motor the relative amount of steps. This function does not enable or disable the motor automatically.

Parameters `steps` – Number of steps to turn the motor.

```
class ILS2TConfig(rpm_max_init: numbers.Integral = 1500, wait_sec_post_enable: Union[int, float] = 1,
                   wait_sec_max_disable: Union[int, float] = 10, wait_sec_post_cannot_disable: Union[int, float] = 1,
                   wait_sec_post_relative_step: Union[int, float] = 2,
                   wait_sec_post_absolute_position: Union[int, float] = 2)
```

Bases: `object`

Configuration for the ILS2T stepper motor device.

`clean_values()`

`force_value(fieldname, value)`

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define `post_force_value` method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field

- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

rpm_max_init: numbers.Integral = 1500

initial maximum RPM for the motor, can be set up to 3000 RPM. The user is allowed to set a new max RPM at runtime using [ILS2T.set_max_rpm\(\)](#), but the value must never exceed this configuration setting.

wait_sec_max_disable: Union[int, float] = 10

wait_sec_post_absolute_position: Union[int, float] = 2

wait_sec_post_CANNOT_DISABLE: Union[int, float] = 1

wait_sec_post_enable: Union[int, float] = 1

wait_sec_post_relative_step: Union[int, float] = 2

exception ILS2TException

Bases: [hvl_ccb.dev.base.DeviceException](#)

Exception to indicate problems with the SE ILS2T stepper motor.

class ILS2TModbusTcpCommunication(configuration)

Bases: [hvl_ccb.comm.modbus_tcp.ModbusTcpCommunication](#)

Specific implementation of Modbus/TCP for the Schneider Electric ILS2T stepper motor.

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

class ILS2TModbusTcpCommunicationConfig(host: Union[str, ipaddress.IPv4Address,

ipaddress.IPv6Address], unit: int = 255, port: int = 502)

Bases: [hvl_ccb.comm.modbus_tcp.ModbusTcpCommunicationConfig](#)

Configuration dataclass for Modbus/TCP communication specific for the Schneider Electric ILS2T stepper motor.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field

- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

unit: int = 255

The unit has to be 255 such that IO scanning mode works.

class ILS2TRegAddr(value)

Bases: enum.IntEnum

Modbus Register Adresses for Schneider Electric ILS2T stepper drive.

ACCESS_ENABLE = 282

FLT_INFO = 15362

FLT_MEM_DEL = 15112

FLT_MEM_RESET = 15114

IO_SCANNING = 6922

JOGN_FAST = 10506

JOGN_SLOW = 10504

POSITION = 7706

RAMP_ACC = 1556

RAMP_DECEL = 1558

RAMP_N_MAX = 1554

RAMP_TYPE = 1574

SCALE = 1550

TEMP = 7200

VOLT = 7198

class ILS2TRegDatatype(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)

Bases: aenum.Enum

Modbus Register Datatypes for Schneider Electric ILS2T stepper drive.

From the manual of the drive:

datatype	byte	min	max
INT8	1 Byte	-128	127
UINT8	1 Byte	0	255
INT16	2 Byte	-32_768	32_767
UINT16	2 Byte	0	65_535
INT32	4 Byte	-2_147_483_648	2_147_483_647
UINT32	4 Byte	0	4_294_967_295
BITS	just 32bits	N/A	N/A

INT32 = (-2147483648, 2147483647)

is_in_range(value: int) → bool

exception IoScanningModeValueError

Bases: *hvl_ccb.dev.se_ilst2t.ILS2TException*


Exception to indicate that the selected IO scanning mode is invalid.

exception ScalingFactorValueError

Bases: *hvl_ccb.dev.se_ilst2t.ILS2TException*

Exception to indicate that a scaling factor value is invalid.

hvl_ccb.dev.sst_luminox

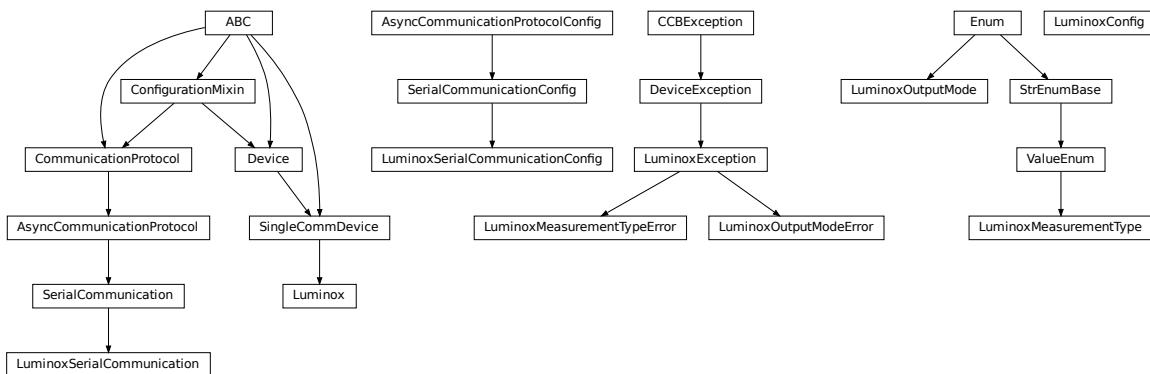

Device class for a SST Luminox Oxygen sensor. This device can measure the oxygen concentration between 0 % and 25 %.

Furthermore, it measures the barometric pressure and internal temperature. The device supports two operating modes: in streaming mode the device measures all parameters every second, in polling mode the device measures only after a query.

Technical specification and documentation for the device can be found at the manufacturer's page: <https://www.sstsensing.com/product/luminox-optical-oxygen-sensors-2/>

class Luminox(com, dev_config=None)

Bases: *hvl_ccb.dev.base.SingleCommDevice*

Luminox oxygen sensor device class.

activate_output(mode: hvl_ccb.dev.sst_luminox.LuminoxOutputMode) → None

activate the selected output mode of the Luminox Sensor. :param mode: polling or streaming

static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

static default_com_cls()

Get the class for the default communication protocol used with this device.

Returns the type of the standard communication protocol for this device

query_polling(measurement: Union[str, hvl_ccb.dev.sst_luminox.LuminoxMeasurementType]) → Union[Dict[Union[str, hvl_ccb.dev.sst_luminox.LuminoxMeasurementType], Union[float, int, str]], float, int, str]

Query a value or values of Luminox measurements in the polling mode, according to a given measurement type.

Parameters **measurement** – type of measurement

Returns value of requested measurement

Raises

- **ValueError** – when a wrong key for LuminoxMeasurementType is provided
- **LuminoxOutputModeError** – when polling mode is not activated
- **LuminoxMeasurementTypeError** – when expected measurement value is not read

read_streaming() → Dict[Union[str, hvl_ccb.dev.sst_luminox.LuminoxMeasurementType], Union[float, int, str]]

Read values of Luminox in the streaming mode. Convert the single string into separate values.

Returns dictionary with *LuminoxMeasurementType.all_measurements_types()* keys and accordingly type-parsed values.

Raises

- **LuminoxOutputModeError** – when streaming mode is not activated
- **LuminoxMeasurementTypeError** – when any of expected measurement values is not read

start() → None

Start this device. Opens the communication protocol.

stop() → None

Stop the device. Closes also the communication protocol.

class LuminoxConfig(wait_sec_post_activate: Union[int, float] = 0.5, wait_sec_trials_activate: Union[int, float] = 0.1, nr_trials_activate: int = 5)

Bases: object

Configuration for the SST Luminox oxygen sensor.

clean_values()**force_value(fieldname, value)**

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field

- **value** – value to assign

is_configdataclass = True

classmethod keys() → Sequence[str]
Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

nr_trials_activate: int = 5

classmethod optional_defaults() → Dict[str, object]
Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]
Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

wait_sec_post_activate: Union[int, float] = 0.5

wait_sec_trials_activate: Union[int, float] = 0.1

exception LuminoxException
Bases: *hvl_ccb.dev.base.DeviceException*
General Exception for Luminox Device.

class LuminoxMeasurementType(*value=<no_arg>*, *names=None*, *module=None*, *type=None*, *start=1*,
boundary=None)
Bases: *hvl_ccb.utils.enum.ValueEnum*
Measurement types for *LuminoxOutputMode.polling*.

The *all_measurements* type will read values for the actual measurement types as given in *LuminoxOutputMode.all_measurements_types()*; it parses multiple single values using regexp's for other measurement types, therefore, no regexp is defined for this measurement type.

all_measurements = 'A'

classmethod all_measurements_types() → Tuple[*hvl_ccb.dev.sst_luminox.LuminoxMeasurementType*, ...]
A tuple of *LuminoxMeasurementType* enum instances which are actual measurements, i.e. not date of manufacture or software revision.

barometric_pressure = 'P'

property command: str

date_of_manufacture = '# 0'

parse_read_measurement_value(*read_txt: str*) → Union[Dict[Union[str, *hvl_ccb.dev.sst_luminox.LuminoxMeasurementType*], Union[float, int, str]], float, int, str]

partial_pressure_o2 = '0'

percent_o2 = '%'

sensor_status = 'e'

serial_number = '# 1'

```

software_revision = '# 2'
temperature_sensor = 'T'

LuminoxMeasurementTypeDict
A typing hint for a dictionary holding LuminoxMeasurementType values. Keys are allowed as strings because LuminoxMeasurementType is of a StrEnumBase type.

alias of Dict[Union[str, LuminoxMeasurementType], Union[float, int, str]]

exception LuminoxMeasurementTypeError
Bases: hvl_ccb.dev.sst_luminox.LuminoxException

Wrong measurement type for requested data

LuminoxMeasurementTypeValue
A typing hint for all possible LuminoxMeasurementType values as read in either streaming mode or in a polling mode with LuminoxMeasurementType.all_measurements.

Beware: has to be manually kept in sync with LuminoxMeasurementType instances cast_type attribute values.

alias of Union[float, int, str]

class LuminoxOutputMode(value)
Bases: enum.Enum

output mode.

polling = 1
streaming = 0

exception LuminoxOutputModeError
Bases: hvl_ccb.dev.sst_luminox.LuminoxException

Wrong output mode for requested data

class LuminoxSerialCommunication(configuration)
Bases: hvl_ccb.comm.serial.SerialCommunication

Specific communication protocol implementation for the SST Luminox oxygen sensor. Already predefines device-specific protocol parameters in config.

static config_cls()
Return the default configdataclass class.

Returns a reference to the default configdataclass class

class LuminoxSerialCommunicationConfig(terminator: bytes = b'\r\n', encoding: str = 'utf-8', encoding_error_handling: str = 'strict', wait_sec_read_text_nonempty: Union[int, float] = 0.5, default_n_attempts_read_text_nonempty: int = 10, port: Union[str, NoneType] = None, baudrate: int = 9600, parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = <SerialCommunicationParity.NONE: 'N'>, stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = <SerialCommunicationStopbits.ONE: 1>, bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = <SerialCommunicationBytesize.EIGHTBITS: 8>, timeout: Union[int, float] = 3)
Bases: hvl_ccb.comm.serial.SerialCommunicationConfig

baudrate: int = 9600
Baudrate for SST Luminox is 9600 baud

```

```
bytesize: Union[int, hvl_ccb.comm.serial.SerialCommunicationBytesize] = 8
    One byte is eight bits long

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define post_force_value method with same signature as this method to do extra processing
after value has been forced on fieldname.
```

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

parity: Union[str, hvl_ccb.comm.serial.SerialCommunicationParity] = 'N'
 SST Luminox does not use parity

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

stopbits: Union[int, hvl_ccb.comm.serial.SerialCommunicationStopbits] = 1
 SST Luminox does use one stop bit

terminator: bytes = b'\r\n'
 The terminator is CR LF

timeout: Union[int, float] = 3
 use 3 seconds timeout as default

hvl_ccb.dev.utils

Poller

```
class Poller(spoll_handler: Callable, polling_delay_sec: Union[int, float] = 0, polling_interval_sec: Union[int,
    float] = 1, polling_timeout_sec: Optional[Union[int, float]] = None)
Bases: object
```

Poller class wrapping *concurrent.futures.ThreadPoolExecutor* which enables passing of results and errors out of the polling thread.

is_polling() → bool

Check if device status is being polled.

Returns *True* when polling thread is set and alive

start_polling() → bool

Start polling.

Returns *True* if was not polling before, *False* otherwise

stop_polling() → bool

Stop polling.

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

Returns *True* if was polling before, *False* otherwise, and last result of the polling function call.


Raises polling function exceptions

wait_for_polling_result()

Wait for until polling function returns a result as well as any exception that might have been raised within a thread.

Returns polling function result

Raises polling function errors

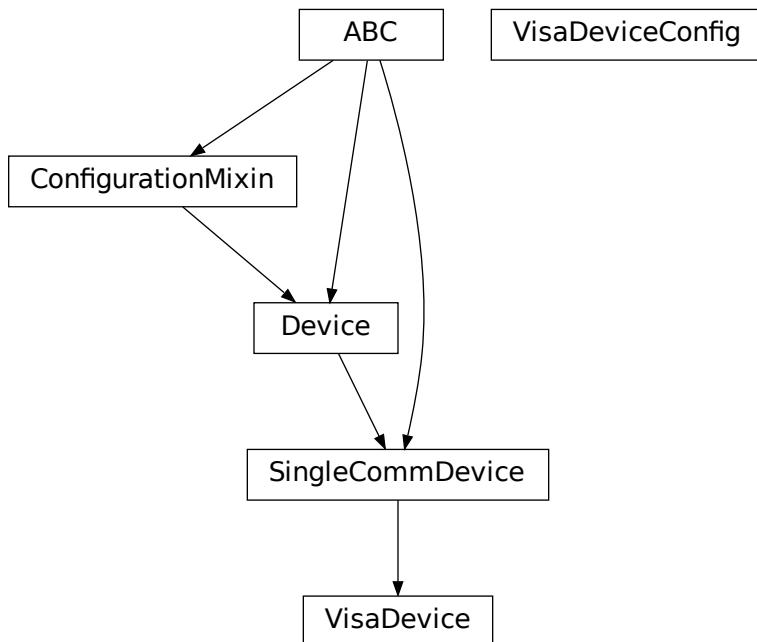hvl_ccb.dev.visa

```
class VisaDevice(com: Union[hvl_ccb.comm.visa.VisaCommunication,
hvl_ccb.comm.visa.VisaCommunicationConfig, dict], dev_config:
Optional[Union[hvl_ccb.dev.visa.VisaDeviceConfig, dict]] = None)
Bases: hvl_ccb.dev.base.SingleCommDevice

Device communicating over the VISA protocol using VisaCommunication.

static config_cls()
    Return the default configdataclass class.

    Returns a reference to the default configdataclass class

static default_com_cls() → Type[hvl_ccb.comm.visa.VisaCommunication]
    Return the default communication protocol for this device type, which is VisaCommunication.

    Returns the VisaCommunication class

get_error_queue() → str
    Read out error queue and logs the error.

    Returns Error string

get_identification() → str
    Queries “*IDN?” and returns the identification string of the connected device.

    Returns the identification string of the connected device

reset() → None
    Send “*RST” and “*CLS” to the device. Typically sets a defined state.

spoll_handler()
    Reads the status byte and decodes it. The status byte STB is defined in IEEE 488.2. It provides a rough
    overview of the instrument status.

    Returns

start() → None
    Start the VisaDevice. Sets up the status poller and starts it.

    Returns

stop() → None
    Stop the VisaDevice. Stops the polling thread and closes the communication protocol.

    Returns

wait_operation_complete(timeout: Optional[float] = None) → bool
    Waits for a operation complete event. Returns after timeout [s] has expired or the operation complete event
    has been caught.

    Parameters timeout – Time in seconds to wait for the event; None for no timeout.

    Returns True, if OPC event is caught, False if timeout expired

class VisaDeviceConfig(spoll_interval: Union[int, float] = 0.5, spoll_start_delay: Union[int, float] = 2)
Bases: hvl_ccb.dev.visa._VisaDeviceConfigDefaultsBase, hvl_ccb.dev.visa.
_VisaDeviceConfigBase

Configdataclass for a VISA device.

force_value(fieldname, value)
    Forces a value to a dataclass field despite the class being frozen.

    NOTE: you can define post_force_value method with same signature as this method to do extra processing
    after value has been forced on fieldname.
```

Parameters

- **fieldname** – name of the field
- **value** – value to assign

classmethod keys() → Sequence[str]

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.


Module contents

Devices subpackage.

hvl_ccb.utils

Submodules

hvl_ccb.utils.conversion_sensor

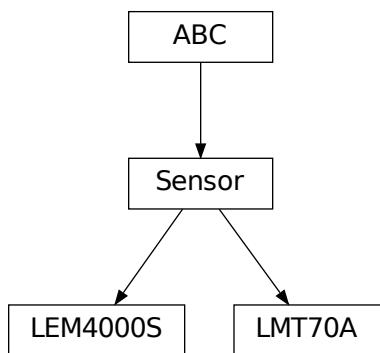

Sensors that are used by the devices implemented in the CCB

class LEM4000S(shunt: float = 1.2, calibration_factor: float = 1)

Bases: [hvl_ccb.utils.conversion_sensor.Sensor](#)

Converts the output voltage (V) to the measured current (A) when using a LEM Current transducer LT 4000-S

```
CONVERSION: ClassVar[int] = 5000
calibration_factor: float = 1
convert(value, **kwargs)
shunt: float = 1.2

class LMT70A(temperature_unit: hvl_ccb.utils.conversion_unit.Temperature = Temperature.C)
Bases: hvl_ccb.utils.conversion_sensor.Sensor

Converts the output voltage (V) to the measured temperature (default °C) when using a TI Precision Analog
Temperature Sensor LMT70(A)

LUT: ClassVar[numpy.ndarray[Any,
numpy.dtype[numpy.typing._generic_alias.ScalarType]]] = array([[-55. , 1.375219],
[-50. , 1.350441], [-40. , 1.300593], [-30. , 1.250398], [-20. , 1.199884], [-10. ,
1.14907], [ 0. , 1.097987], [ 10. , 1.046647], [ 20. , 0.99505], [ 30. ,
0.943227], [ 40. , 0.891178], [ 50. , 0.838882], [ 60. , 0.78636], [ 70. ,
0.733608], [ 80. , 0.680654], [ 90. , 0.62749], [100. , 0.574117], [110. ,
0.520551], [120. , 0.46676], [130. , 0.412739], [140. , 0.358164], [150. ,
0.302785]])]


convert(value, **kwargs)
temperature_unit: hvl_ccb.utils.conversion_unit.Temperature = 'C'

class Sensor
Bases: abc.ABC

The BaseClass ‘Sensor’ is designed as a parent for all Sensors. Each attribute must be added to ‘__setattr__’, so
that the value is verified each time the value is changed. It is important to mark attributes that should be constant
with ‘typing.ClassVar[...]’. Together with ‘super().__setattr__(name, value)’, this guarantees that the values are
protected and cannot be altered by the user.

abstract convert(value, **kwargs)
```

hvl_ccb.utils.conversion_unit

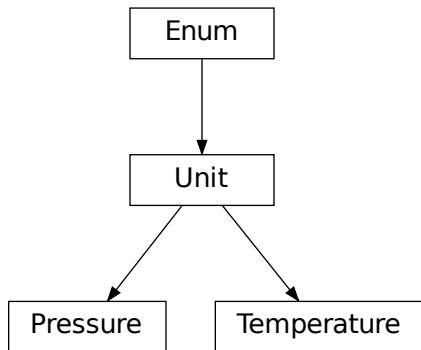

Unit conversion, within in the same group of units, for example Kelvin <-> Celsius

```
class Pressure(value)
Bases: hvl_ccb.utils.conversion_unit.Unit

An enumeration.

ATM = 'atm'
ATMOSPHERE = 'atm'
BAR = 'bar'
MILLIMETER_MERCURY = 'mmHg'
MMHG = 'mmHg'
PA = 'Pa'
PASCAL = 'Pa'
POUNDS_PER_SQUARE_INCH = 'psi'
PSI = 'psi'
TORR = 'torr'

class Temperature(value)
Bases: hvl_ccb.utils.conversion_unit.Unit

An enumeration.

C = 'C'
CELSIUS = 'C'
F = 'F'
FAHRENHEIT = 'F'
K = 'K'
KELVIN = 'K'


class Unit(value)
Bases: enum.Enum

An enumeration.

abstract classmethod convert(value, **kwargs)
```

preserve_type(func)

This wrapper preserves the first order type of the input. Upto now the type of the data stored in a list, tuple, array or dict is not preserved. Integer will be converted to float!

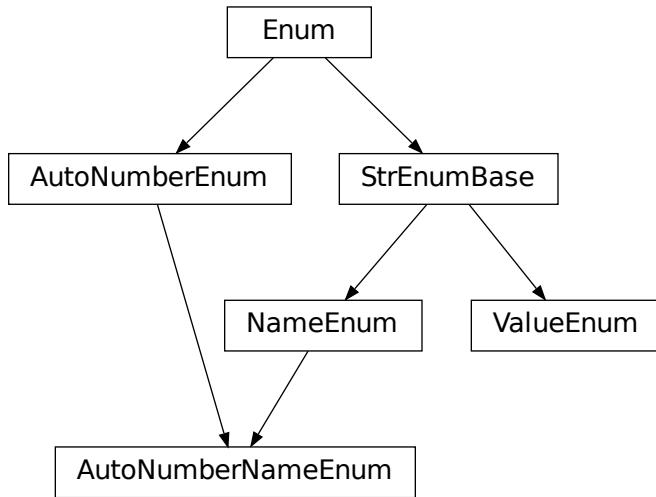hvl_ccb.utils.enum

```
class AutoNumberNameEnum(value=<no_arg>, names=None, module=None, type=None, start=1,  
boundary=None)
```

Bases: [hvl_ccb.utils.enum.NameEnum](#), [aenum.AutoNumberEnum](#)

Auto-numbered enum with names used as string representation, and with lookup and equality based on this representation.

```
class NameEnum(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: hvl\_ccb.utils.enum.StrEnumBase
```

Enum with names used as string representation, and with lookup and equality based on this representation.

```
class StrEnumBase(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: aenum.Enum
```

String representation-based equality and lookup.

```
class ValueEnum(value=<no_arg>, names=None, module=None, type=None, start=1, boundary=None)  
Bases: hvl\_ccb.utils.enum.StrEnumBase
```

Enum with string representation of values used as string representation, and with lookup and equality based on this representation.

Attention: to avoid errors, best use together with *unique* enum decorator.

hvl_ccb.utils.typing

Additional Python typing module utilities

ConvertibleTypes

Typing hint for data type that can be used in conversion

alias of Union[int, float, List[Union[int, float]], Tuple[Union[int, float]], Dict[str, Union[int, float]], numpy.ndarray]

Number

Typing hint auxiliary for a Python base number types: *int* or *float*.

alias of Union[int, float]

check_generic_type(value, type_, name='instance')

Check if *value* is of a generic type *type_*. Raises *TypeError* if it's not.

Parameters

- **name** – name to report in case of an error
- **value** – value to check
- **type** – generic type to check against

is_generic_type_hint(type_)

Check if class is a generic type, for example *Union[int, float]*, *List[int]*, or *List*.

Parameters **type** – type to check

hvl_ccb.utils.validation

validate_and_resolve_host(host: Union[str, ipaddress.IPv4Address, ipaddress.IPv6Address], logger: Optional[logging.Logger] = None) → str

validate_bool(x_name: str, x: object, logger: Optional[logging.Logger] = None) → None

Validate if given input *x* is a *bool*.

Parameters

- **x_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as boolean
- **logger** – logger of the calling submodule

Raises **TypeError** – when the validated input does not have boolean type

validate_number(x_name: str, x: object, limits: Optional[Tuple] = (None, None), number_type: Union[Type[Union[int, float]], Tuple[Type[Union[int, float]], ...]] = (<class 'int'>, <class 'float'>), logger: Optional[logging.Logger] = None) → None

Validate if given input *x* is a number of given *number_type* type, with value between given *limits[0]* and *limits[1]* (inclusive), if not *None*. For array-like objects (npt.NDArray, List, Tuple, Dict) it is checked if all elements are within the limits and have the correct type.

Parameters

- **x_name** – string name of the validate input, use for the error message
- **x** – an input object to validate as number of given type within given range
- **logger** – logger of the calling submodule

- **limits** – [lower, upper] limit, with *None* denoting no limit: $[-\infty, +\infty]$
- **number_type** – expected type or tuple of types of a number, by default (*int*, *float*)

Raises


- **TypeError** – when the validated input does not have expected type
- **ValueError** – when the validated input has correct number type but is not within given range

validate_tcp_port(*port*: *int*, *logger*: *Optional[logging.Logger]* = *None*)

Module contents

4.1.2 Submodules

[hvl_ccb.configuration](#)

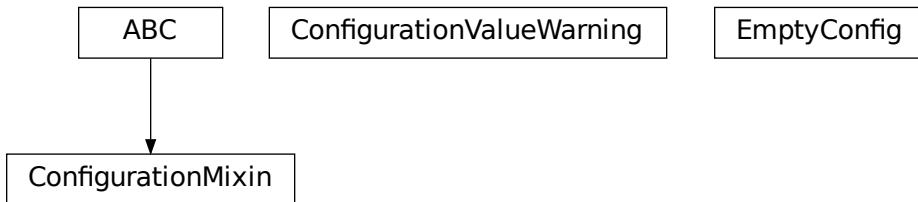

providing classes for handling configuration for communication protocols and devices.

Facilities

class ConfigurationMixin(*configuration*)

Bases: abc.ABC

Mixin providing configuration to a class.

property config

ConfigDataclass property.

Returns the configuration

abstract static config_cls()

Return the default configdataclass class.

Returns a reference to the default configdataclass class

configuration_save_json(*path*: str) → None

Save current configuration as JSON file.

Parameters *path* – path to the JSON file.

classmethod from_json(*filename*: str)

Instantiate communication protocol using configuration from a JSON file.

Parameters *filename* – Path and filename to the JSON configuration

exception ConfigurationValueWarning

Bases: UserWarning

User warnings category for values of @configdataclass fields.

class EmptyConfig

Bases: object

Empty configuration dataclass.

clean_values()

Cleans and enforces configuration values. Does nothing by default, but may be overridden to add custom configuration value checks.

force_value(fieldname, value)

Forces a value to a dataclass field despite the class being frozen.

NOTE: you can define *post_force_value* method with same signature as this method to do extra processing after *value* has been forced on *fieldname*.

Parameters

- **fieldname** – name of the field
- **value** – value to assign

is_configdataclass = True**classmethod keys() → Sequence[str]**

Returns a list of all configdataclass fields key-names.

Returns a list of strings containing all keys.

classmethod optional_defaults() → Dict[str, object]

Returns a list of all configdataclass fields, that have a default value assigned and may be optionally specified on instantiation.

Returns a list of strings containing all optional keys.

classmethod required_keys() → Sequence[str]

Returns a list of all configdataclass fields, that have no default value assigned and need to be specified on instantiation.

Returns a list of strings containing all required keys.

configdataclass(direct_decoration=None, frozen=True)

Decorator to make a class a configdataclass. Types in these dataclasses are enforced. Implement a function clean_values(self) to do additional checking on value ranges etc.

It is possible to inherit from a configdataclass and re-decorate it with @configdataclass. In a subclass, default values can be added to existing fields. Note: adding additional non-default fields is prone to errors, since the order has to be respected through the whole chain (first non-default fields, only then default-fields).

Parameters **frozen** – defaults to True. False allows to later change configuration values. Attention:
if configdataclass is not frozen and a value is changed, typing is not enforced anymore!

[hvl_ccb.exception](#)


```
CCBException
```

Introduce a common base exception for the CCB

```
exception CCBException
```

Bases: Exception

[hvl_ccb.experiment_manager](#)

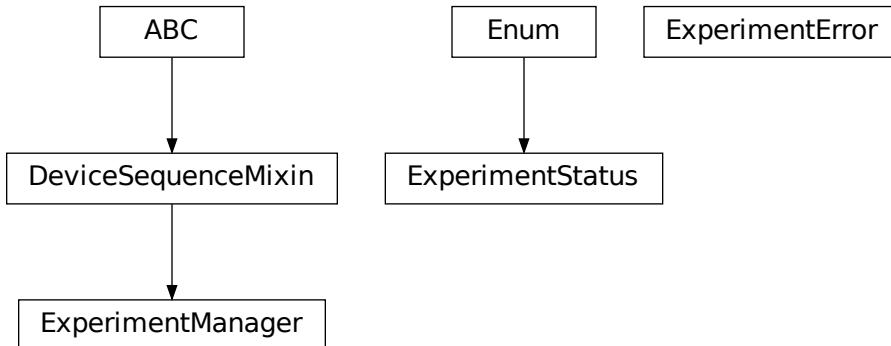

Main module containing the top level ExperimentManager class. Inherit from this class to implement your own experiment functionality in another project and it will help you start, stop and manage your devices.

```
exception ExperimentError
```

Bases: Exception

Exception to indicate that the current status of the experiment manager is on ERROR and thus no operations can be made until reset.

```
class ExperimentManager(*args, **kwargs)
```

Bases: [hvl_ccb.dev.base.DeviceSequenceMixin](#)

Experiment Manager can start and stop communication protocols and devices. It provides methods to queue commands to devices and collect results.

```
add_device(name: str, device: hvl_ccb.dev.base.Device) → None
```

Add a new device to the manager. If the experiment is running, automatically start the device. If a device with this name already exists, raise an exception.

Parameters

- **name** – is the name of the device.

- **device** – is the instantiated Device object.

Raises *DeviceExistingException* –

devices_failed_start: `Dict[str, hvl_ccb.dev.base.Device]`

Dictionary of named device instances from the sequence for which the most recent *start()* attempt failed.

Empty if *stop()* was called last; cf. *devices_failed_stop*.

devices_failed_stop: `Dict[str, hvl_ccb.dev.base.Device]`

Dictionary of named device instances from the sequence for which the most recent *stop()* attempt failed.

Empty if *start()* was called last; cf. *devices_failed_start*.

finish() → None

Stop experimental setup, stop all devices.

is_error() → bool

Returns true, if the status of the experiment manager is *error*.

Returns True if on error, false otherwise

is_finished() → bool

Returns true, if the status of the experiment manager is *finished*.

Returns True if finished, false otherwise

is_running() → bool

Returns true, if the status of the experiment manager is *running*.

Returns True if running, false otherwise

run() → None

Start experimental setup, start all devices.

start() → None

Alias for ExperimentManager.run()

property status: `hvl_ccb.experiment_manager.ExperimentStatus`

Get experiment status.

Returns experiment status enum code.

stop() → None

Alias for ExperimentManager.finish()

class ExperimentStatus(*value*)

Bases: `enum.Enum`

Enumeration for the experiment status

ERROR = 5

FINISHED = 4

FINISHING = 3

INITIALIZED = 0

INITIALIZING = -1

RUNNING = 2

STARTING = 1

4.1.3 Module contents

Top-level package for HVL Common Code Base.

CONTRIBUTING

Contributions are welcome, and they are greatly appreciated! Every little bit helps, and credit will always be given.

You can contribute in many ways:

5.1 Types of Contributions

5.1.1 Report Bugs

Report bugs at https://gitlab.com/ethz_hvl/hvl_ccb/issues.

If you are reporting a bug, please include:

- Your operating system name and version.
- Any details about your local setup that might be helpful in troubleshooting.
- Detailed steps to reproduce the bug.

5.1.2 Fix Bugs

Look through the GitLab issues for bugs. Anything tagged with “bug” and “help wanted” is open to whoever wants to implement it.

5.1.3 Implement Features

Look through the GitLab issues for features. Anything tagged with “enhancement” and “help wanted” is open to whoever wants to implement it.

5.1.4 Write Documentation

HVL Common Code Base could always use more documentation, whether as part of the official HVL Common Code Base docs, in docstrings, or even on the web in blog posts, articles, and such.

5.1.5 Submit Feedback

The best way to send feedback is to file an issue at https://gitlab.com/ethz_hvl/hvl_ccb/issues.

If you are proposing a feature:

- Explain in detail how it would work.
- Keep the scope as narrow as possible, to make it easier to implement.
- Remember that this is a volunteer-driven project, and that contributions are welcome :)

5.2 Get Started!

Ready to contribute? Here's how to set up *hvl_ccb* for local development.

1. Clone *hvl_ccb* repo from GitLab.

```
$ git clone git@gitlab.com:ethz_hvl/hvl_ccb.git
```

2. Install your local copy into a virtualenv. Assuming you have `virtualenvwrapper` installed, this is how you set up your fork for local development:

```
$ mkvirtualenv hvl_ccb
$ cd hvl_ccb/
$ pip install -e .[all]
$ pip install -r requirements_dev.txt
```

3. Create a branch for local development:

```
$ git checkout -b name-of-your-bugfix-or-feature
```

Now you can make your changes locally.

4. When you're done making changes, check that your changes pass flake8 and the tests, including testing other Python versions with tox:

```
$ flake8 hvl_ccb tests
$ python setup.py test or py.test
$ tox
```

To get flake8 and tox, just pip install them into your virtualenv. You can also use the provided make-like shell script to run flake8 and tests:

```
$ ./make.sh lint
$ ./make.sh test
```

5. Commit your changes and push your branch to GitLab:

```
$ git add .
$ git commit -m "Your detailed description of your changes."
$ git push origin name-of-your-bugfix-or-feature
```

6. Submit a merge request through the GitLab website.

5.3 Merge Request Guidelines

Before you submit a merge request, check that it meets these guidelines:

1. The merge request should include tests.
2. If the merge request adds functionality, the docs should be updated. Put your new functionality into a function with a docstring, and add the feature to the list in README.rst.
3. The merge request should work for Python 3.7. Check https://gitlab.com/ethz_hvl/hvl_ccb/merge_requests and make sure that the tests pass for all supported Python versions.

5.4 Tips

- To run tests from a single file:

```
$ py.test tests/test_hvl_ccb.py
```

or a single test function:

```
$ py.test tests/test_hvl_ccb.py::test_command_line_interface
```

- If your tests are slow, profile them using the pytest-profiling plugin:

```
$ py.test tests/test_hvl_ccb.py --profile
```

or for a graphical overview (you need a SVG image viewer):

```
$ py.test tests/test_hvl_ccb.py --profile-svg  
$ open prof/combined.svg
```

- To add dependency, edit appropriate *requirements variable in the setup.py file and re-run:

```
$ python setup.py develop
```

- To generate a PDF version of the Sphinx documentation instead of HTML use:

```
$ rm -rf docs/hvl_ccb.rst docs/modules.rst docs/_build && sphinx-apidoc -o docs/hvl_ccb && python -msphinx -M latexpdf docs/ docs/_build
```

This command can also be run through the make-like shell script:

```
$ ./make.sh docs-pdf
```

This requires a local installation of a LaTeX distribution, e.g. MikTeX.

5.5 Deploying

A reminder for the maintainers on how to deploy. Create `release-N.M.K` branch. Make sure all your changes are committed. Update or create entry in `HISTORY.rst` file, and, if applicable, update `AUTHORS.rst` file, update features tables in `README.rst` file, and update API docs:

```
$ make docs
```

Commit all of the above, except for the `docs/hvl_ccb.dev.picotech_pt104.rst`, and then run:

```
$ bumpversion patch # possible: major / minor / patch  
$ git push  
$ git push --tags
```

Go to <https://readthedocs.org/projects/hvlccb/builds/> and check if RTD docs build for the pushed tag passed.

Wait for the CI pipeline to finish successfully; afterwards, run a release check:

```
$ make release-check
```

Finally, prepare and push a release:

```
$ make release
```

Merge the release branch into master and devel branches with `--no-ff` flag and delete the release branch:

```
$ git switch master  
$ git merge --no-ff release-N.M.K  
$ git push  
$ git switch devel  
$ git merge --no-ff release-N.M.K  
$ git push  
$ git push --delete origin release-N.M.K  
$ git branch --delete release-N.M.K
```

Finally, prepare GitLab release and cleanup the corresponding milestone:

1. go to https://gitlab.com/ethz_hvl/hvl_ccb/-/tags/, select the latest release tag, press “Edit release notes” and add the release notes (copy a corresponding entry from `HISTORY.rst` file with formatting adjusted from ReStructuredText to Markdown); press “Save changes”;
2. go to https://gitlab.com/ethz_hvl/hvl_ccb/-/releases, select the latest release, press “Edit this release” and under “Milestones” select the corresponding milestone; press “Save changes”;
3. go to https://gitlab.com/ethz_hvl/hvl_ccb/-/milestones, make sure that it is 100% complete (otherwise, create a next patch-level milestone and assign it to the ongoing Issues and Merge Requests therein); press “Close Milestone”.

CREDITS

6.1 Maintainers

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>

6.2 Authors

- Mikołaj Rybiński <mikolaj.rybinski@id.ethz.ch>
- David Gruber <dev@davidgruber.ch>
- Henrik Menne <henrik.menne@eeh.ee.ethz.ch>
- Alise Chachereau <chachereau@eeh.ee.ethz.ch>
- Henning Janssen <janssen@eeh.ee.ethz.ch>
- David Taylor <dtaylor@ethz.ch>
- Joseph Engelbrecht <engelbrecht@eeh.ee.ethz.ch>

6.3 Contributors

- Luca Nembrini <lucane@student.ethz.ch>
- Maria Del <maria.del@id.ethz.ch>
- Raphael Faerber <raphael.faerber@eeh.ee.ethz.ch>
- Ruben Stadler <rstadler@student.ethz.ch>
- Hanut Vemulapalli <vemulapalli@eeh.ee.ethz.ch>

HISTORY

7.1 0.9.0 (2022-01-07)

- New device: Highland T560 digital delay and pulse generator over Telnet.
- **Rework of the Technix Capacitor Charger.**
 - Moved into a separate sub-package
 - NEW import over `import hvl_ccb.dev.technix as XXX`
 - Slightly adapted behaviour
- Add `validate_tcp_port` to validate port number.
- **Add `validate_and_resolve_host` to validate and resolve host names and IPs.**
 - Remove requirement IPy
- Add a unified CCB Exception schema for all devices and communication protocols.
- Add data conversion functions to README.
- Update CI and devel images from Debian 10 buster to Debian 11 bullseye.
- Fix typing due to numpy update.
- Fix incorrect overloading of `clean_values()` in classes of type `XCommunicationConfig`.

7.2 0.8.5 (2021-11-05)

- Added arbitrary waveform for TiePie signal generation, configurable via `dev.tiepie.generator.TiePieGeneratorConfig.waveform` property.
- In `utils.conversion_sensor`: improvements for class constants; removed SciPy dependency.
- Added Python 3.10 support.

7.3 0.8.4 (2021-10-22)

- `utils.validation.validate_number` extension to handle NumPy arrays and array-like objects.
- `utils.conversion_unit` utility classes handle correctly `NamedTuple` instances.
- `utils.conversion_sensor` and `utils.conversion_unit` code simplification (no `transfer_function_order` attribute) and cleanups.
- Fixed incorrect error logging in `configuration.configdataclass`.
- `comm.telnet.TelnetCommunication` tests fixes for local run errors.

7.4 0.8.3 (2021-09-27)

- New data conversion functions in `utils.conversion_sensor` and `utils.conversion_unit` modules. Note: to use these functions you must install `hvl_ccb` with extra requirement, either `hvl_ccb[conversion]` or `hvl_ccb[all]`.
- Improved documentation with respect to installation of external libraries.

7.5 0.8.2 (2021-08-27)

- New functionality in `dev.labjack.LabJack`:
 - configure clock and send timed pulse sequences
 - set DAC/analog output voltage
- Bugfix: ignore random bits sent by to `dev.newport.NewportSMC100PP` controller during start-up/powering-up.

7.6 0.8.1 (2021-08-13)

- Add Python version check (min version error; max version warning).
- Daily checks for upstream dependencies compatibility and devel environment improvements.

7.7 0.8.0 (2021-07-02)

- TCP communication protocol.
- Lauda PRO RP 245 E circulation thermostat device over TCP.
- Pico Technology PT-104 Platinum Resistance Data Logger device as a wrapper of the Python bindings for the PicoSDK.
- In `com.visa.VisaCommunication`: periodic status polling when VISA/TCP keep alive connection is not supported by a host.

7.8 0.7.1 (2021-06-04)

- New `utils.validation` submodule with `validate_bool` and `validate_number` utilities extracted from internal use within a `dev.tiepie` subpackage.
- In `comm.serial.SerialCommunication`:
 - strict encoding errors handling strategy for subclasses,
 - user warning for a low communication timeout value.

7.9 0.7.0 (2021-05-25)

- The `dev.tiepie` module was splitted into a subpackage with, in particular, submodules for each of the device types – oscilloscope, generator, and i2c – and with backward-incompatible direct imports from the submodules.
- In `dev.technix`:
 - fixed communication crash on nested status byte query;
 - added enums for GET and SET register commands.
- Further minor logging improvements: added missing module level logger and removed some error logs in `except` blocks used for a flow control.
- In `examples/` folder renamed consistently all the examples.
- In API documentation: fix incorrect links mapping on inheritance diagrams.

7.10 0.6.1 (2021-05-08)

- In `dev.tiepie`:
 - dynamically set oscilloscope's channel limits in `OscilloscopeChannelParameterLimits`: `input_range` and `trigger_level_abs`, incl. update of latter on each change of `input_range` value of a `TiePieOscilloscopeChannelConfig` instances;
 - quick fix for opening of combined instruments by disabling `OscilloscopeParameterLimits`. `trigger_delay` (an advanced feature);
 - enable automatic devices detection to be able to find network devices with `TiePieOscilloscope`. `list_devices()`.
- Fix `examples/example_labjack.py`.
- Improved logging: consistently use module level loggers, and always log exception tracebacks.
- Improve API documentation: separate pages per modules, each with an inheritance diagram as an overview.

7.11 0.6.0 (2021-04-23)

- Technix capacitor charger using either serial connection or Telnet protocol.
- **Extensions, improvements and fixes in existing devices:**
 - In `dev.tiepie.TiePieOscilloscope`:
 - * redesigned measurement start and data collection API, incl. time out argument, with no/infinite time out option;
 - * trigger allows now a no/infinite time out;
 - * record length and trigger level were fixed to accept, respectively, floating point and integer numbers;
 - * fixed resolution validation bug;
 - `dev.heinzinger.HeinzingerDI` and `dev.rs_rto1024.RTO1024` instances are now resilient to multiple `stop()` calls.
 - In `dev.crylas.CryLasLaser`: default configuration timeout and polling period were adjusted;
 - Fixed PSI9080 example script.
- **Package and source code improvements:**
 - Update to backward-incompatible `pyvisa-py>=0.5.2`. Developers, do update your local development environments!
 - External libraries, like LibTiePie SDK or LJM Library, are now not installed by default; they are now extra installation options.
 - Added Python 3.9 support.
 - Improved number formatting in logs.
 - Typing improvements and fixes for `mypy>=0.800`.

7.12 0.5.0 (2020-11-11)

- TiePie USB oscilloscope, generator and I2C host devices, as a wrapper of the Python bindings for the LibTiePie SDK.
- a FuG Elektronik Power Supply (e.g. Capacitor Charger HCK) using the built-in ADDAT controller with the Probus V protocol over a serial connection
- All devices polling status or measurements use now a `dev.utils.Poller` utility class.
- **Extensions and improvements in existing devices:**
 - In `dev.rs_rto1024.RTO1024`: added Channel state, scale, range, position and offset accessors, and measurements activation and read methods.
 - In `dev.sst_luminox.Luminox`: added querying for all measurements in polling mode, and made output mode activation more robust.
 - In `dev.newport.NewportSMC100PP`: an error-prone `wait_until_move_finished` method of replaced by a fixed waiting time, device operations are now robust to a power supply cut, and device restart is not required to apply a start configuration.
- **Other minor improvements:**

- Single failure-safe starting and stopping of devices sequenced via `dev.base.DeviceSequenceMixin`.
- Moved `read_text_nonempty` up to `comm.serial.SerialCommunication`.
- Added development Dockerfile.
- Updated package and development dependencies: `pymodbus`, `pytest-mock`.

7.13 0.4.0 (2020-07-16)

- **Significantly improved new Supercube device controller:**
 - more robust error-handling,
 - status polling with generic Poller helper,
 - messages and status boards.
 - tested with a physical device,
- Improved OPC UA client wrapper, with better error handling, incl. re-tries on `concurrent.futures.TimeoutError`.
- SST Luminox Oxygen sensor device controller.
- **Backward-incompatible changes:**
 - `CommunicationProtocol.access_lock` has changed type from `threading.Lock` to `threading.RLock`.
 - `ILS2T.relative_step` and `ILS2T.absolute_position` are now called, respectively, `ILS2T.write_relative_step` and `ILS2T.write_absolute_position`.
- **Minor bugfixes and improvements:**
 - fix use of max resolution in `Labjack.set_ain_resolution()`,
 - resolve ILS2T devices relative and absolute position setters race condition,
 - added acoustic horn function in the 2015 Supercube.
- **Toolchain changes:**
 - add Python 3.8 support,
 - drop pytest-runner support,
 - ensure compatibility with `labjack_ljm` 2019 version library.

7.14 0.3.5 (2020-02-18)

- Fix issue with reading integers from LabJack LJM Library (device's product ID, serial number etc.)
- Fix development requirements specification (tox version).

7.15 0.3.4 (2019-12-20)

- New devices using serial connection:
 - Heinzinger Digital Interface I/II and a Heinzinger PNC power supply
 - Q-switched Pulsed Laser and a laser attenuator from CryLas
 - Newport SMC100PP single axis motion controller for 2-phase stepper motors
 - Pfeiffer TPG controller (TPG 25x, TPG 26x and TPG 36x) for Compact pressure Gauges
- PEP 561 compatibility and related corrections for static type checking (now in CI)
- Refactorings:
 - Protected non-thread safe read and write in communication protocols
 - Device sequence mixin: start/stop, add/rm and lookup
 - `.format()` to f-strings
 - more enumerations and a quite some improvements of existing code
- Improved error docstrings (`:raises:` annotations) and extended tests for errors.

7.16 0.3.3 (2019-05-08)

- Use PyPI labjack-ljm (no external dependencies)

7.17 0.3.2 (2019-05-08)

- INSTALLATION.rst with LJMPython prerequisite info

7.18 0.3.1 (2019-05-02)

- readthedocs.org support

7.19 0.3 (2019-05-02)

- Prevent an automatic close of VISA connection when not used.
- Rhode & Schwarz RTO 1024 oscilloscope using VISA interface over `TCP::INSTR`.
- Extended tests incl. messages sent to devices.
- Added Supercube device using an OPC UA client
- Added Supercube 2015 device using an OPC UA client (for interfacing with old system version)

7.20 0.2.1 (2019-04-01)

- Fix issue with LJMPython not being installed automatically with setuptools.

7.21 0.2.0 (2019-03-31)

- LabJack LJM Library communication wrapper and LabJack device.
- Modbus TCP communication protocol.
- Schneider Electric ILS2T stepper motor drive device.
- Elektro-Automatik PSI9000 current source device and VISA communication wrapper.
- Separate configuration classes for communication protocols and devices.
- Simple experiment manager class.

7.22 0.1.0 (2019-02-06)

- Communication protocol base and serial communication implementation.
- Device base and MBW973 implementation.

**CHAPTER
EIGHT**

INDICES AND TABLES

- genindex
- modindex
- search

PYTHON MODULE INDEX

h

hvl_ccb, 194

hvl_ccb.comm, 32

hvl_ccb.comm.base, 11

hvl_ccb.comm.labjack_ljm, 15

hvl_ccb.comm.modbus_tcp, 18

hvl_ccb.comm.opc, 20

hvl_ccb.comm.serial, 22

hvl_ccb.comm.tcp, 26

hvl_ccb.comm.telnet, 27

hvl_ccb.comm.visa, 29

hvl_ccb.configuration, 190

hvl_ccb.dev, 185

hvl_ccb.dev.base, 95

hvl_ccb.dev.crylas, 97

hvl_ccb.dev.ea_psi9000, 107

hvl_ccb.dev.fug, 111

hvl_ccb.dev.heinzinger, 122

hvl_ccb.dev.highland_t560, 37

hvl_ccb.dev.highland_t560.base, 32

hvl_ccb.dev.highland_t560.channel, 34

hvl_ccb.dev.highland_t560.device, 35

hvl_ccb.dev.labjack, 128

hvl_ccb.dev.lauda, 134

hvl_ccb.dev.mbw973, 141

hvl_ccb.dev.newport, 144

hvl_ccb.dev.pfeiffer_tpg, 158

hvl_ccb.dev.rs_rto1024, 164

hvl_ccb.dev.se_ilis2t, 172

hvl_ccb.dev.sst_luminox, 178

hvl_ccb.dev.supercube, 61

hvl_ccb.dev.supercube.base, 37

hvl_ccb.dev.supercube.constants, 43

hvl_ccb.dev.supercube.typ_a, 57

hvl_ccb.dev.supercube.typ_b, 60

hvl_ccb.dev.supercube2015, 76

hvl_ccb.dev.supercube2015.base, 61

hvl_ccb.dev.supercube2015.constants, 67

hvl_ccb.dev.supercube2015.typ_a, 74

hvl_ccb.dev.technix, 82

hvl_ccb.dev.technix.base, 77

hvl_ccb.dev.technix.device, 80

hvl_ccb.dev.tiepie, 94
hvl_ccb.dev.tiepie.base, 83
hvl_ccb.dev.tiepie.channel, 85
hvl_ccb.dev.tiepie.device, 87
hvl_ccb.dev.tiepie.generator, 88
hvl_ccb.dev.tiepie.i2c, 90
hvl_ccb.dev.tiepie.oscilloscope, 91
hvl_ccb.dev.tiepie.utils, 93
hvl_ccb.dev.utils, 182
hvl_ccb.dev.visa, 183
hvl_ccb.exception, 192
hvl_ccb.experiment_manager, 192
hvl_ccb.utils, 190
hvl_ccb.utils.conversion_sensor, 185
hvl_ccb.utils.conversion_unit, 186
hvl_ccb.utils.enum, 188
hvl_ccb.utils.typing, 189
hvl_ccb.utils.validation, 189

INDEX

A

A (*HeinzingerPNC.UnitCurrent attribute*), 125
A (*SupercubeOpcEndpoint attribute*), 57, 74
ABSOLUTE (*TiePieOscilloscopeTriggerLevelMode attribute*), 87
ABSOLUTE_POSITION (*ILS2T.ActionsPtp attribute*), 172
AC (*NewportConfigCommands attribute*), 144
AC_DoubleStage_150kV (*PowerSetup attribute*), 55, 72
AC_DoubleStage_200kV (*PowerSetup attribute*), 55, 72
AC_SingleStage_100kV (*PowerSetup attribute*), 55, 72
AC_SingleStage_50kV (*PowerSetup attribute*), 55, 72
ACA (*TiePieOscilloscopeChannelCoupling attribute*), 86
acceleration (*NewportSMC100PPConfig attribute*), 152
ACCESS_ENABLE (*ILS2TRegAddr attribute*), 177
access_lock (*CommunicationProtocol attribute*), 14
ACTION_JOG_VALUE (*ILS2T attribute*), 172
activate_clock_output() (*T560 method*), 35
activate_measurements() (*RTO1024 method*), 165
activate_output() (*Luminox method*), 178
activated (*BreakdownDetection attribute*), 49, 68
ACTIVE (*CryLasLaser.AnswersStatus attribute*), 100
active (*OpcControl attribute*), 55
ACTIVE_HIGH (*Polarity attribute*), 33
ACTIVE_LOW (*Polarity attribute*), 33
actualsetvalue (*FuGProbusVSetRegisters property*), 119
ACV (*TiePieOscilloscopeChannelCoupling attribute*), 86
adc_mode (*FuGProbusVMonitorRegisters property*), 119
add_device() (*DeviceSequenceMixin method*), 95
add_device() (*ExperimentManager method*), 192
ADDR_INCORRECT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
address (*NewportSMC100PPConfig attribute*), 152
address (*VisaCommunicationConfig property*), 31
address() (*VisaCommunicationConfig.InterfaceType method*), 30
ADMODE (*FuGProbusIVCommands attribute*), 116
Alarm0 (*AlarmText attribute*), 67
Alarm1 (*Alarms attribute*), 44
Alarm1 (*AlarmText attribute*), 43, 67

Alarm10 (*Alarms attribute*), 44
Alarm10 (*AlarmText attribute*), 43, 67
Alarm100 (*Alarms attribute*), 44
Alarm101 (*Alarms attribute*), 45
Alarm102 (*Alarms attribute*), 45
Alarm103 (*Alarms attribute*), 45
Alarm104 (*Alarms attribute*), 45
Alarm105 (*Alarms attribute*), 45
Alarm106 (*Alarms attribute*), 45
Alarm107 (*Alarms attribute*), 45
Alarm108 (*Alarms attribute*), 45
Alarm109 (*Alarms attribute*), 45
Alarm11 (*Alarms attribute*), 45
Alarm11 (*AlarmText attribute*), 43, 67
Alarm110 (*Alarms attribute*), 45
Alarm111 (*Alarms attribute*), 45
Alarm112 (*Alarms attribute*), 45
Alarm113 (*Alarms attribute*), 45
Alarm114 (*Alarms attribute*), 45
Alarm115 (*Alarms attribute*), 45
Alarm116 (*Alarms attribute*), 45
Alarm117 (*Alarms attribute*), 45
Alarm118 (*Alarms attribute*), 45
Alarm119 (*Alarms attribute*), 45
Alarm12 (*Alarms attribute*), 45
Alarm12 (*AlarmText attribute*), 43, 67
Alarm120 (*Alarms attribute*), 45
Alarm121 (*Alarms attribute*), 45
Alarm122 (*Alarms attribute*), 45
Alarm123 (*Alarms attribute*), 45
Alarm124 (*Alarms attribute*), 45
Alarm125 (*Alarms attribute*), 45
Alarm126 (*Alarms attribute*), 45
Alarm127 (*Alarms attribute*), 45
Alarm128 (*Alarms attribute*), 45
Alarm129 (*Alarms attribute*), 45
Alarm13 (*Alarms attribute*), 45
Alarm13 (*AlarmText attribute*), 43, 67
Alarm130 (*Alarms attribute*), 45
Alarm131 (*Alarms attribute*), 45
Alarm132 (*Alarms attribute*), 45
Alarm133 (*Alarms attribute*), 45

Alarm134 (*Alarms attribute*), 46
Alarm135 (*Alarms attribute*), 46
Alarm136 (*Alarms attribute*), 46
Alarm137 (*Alarms attribute*), 46
Alarm138 (*Alarms attribute*), 46
Alarm139 (*Alarms attribute*), 46
Alarm14 (*Alarms attribute*), 46
Alarm14 (*AlarmText attribute*), 43, 67
Alarm140 (*Alarms attribute*), 46
Alarm141 (*Alarms attribute*), 46
Alarm142 (*Alarms attribute*), 46
Alarm143 (*Alarms attribute*), 46
Alarm144 (*Alarms attribute*), 46
Alarm145 (*Alarms attribute*), 46
Alarm146 (*Alarms attribute*), 46
Alarm147 (*Alarms attribute*), 46
Alarm148 (*Alarms attribute*), 46
Alarm149 (*Alarms attribute*), 46
Alarm15 (*Alarms attribute*), 46
Alarm15 (*AlarmText attribute*), 43
Alarm150 (*Alarms attribute*), 46
Alarm151 (*Alarms attribute*), 46
Alarm16 (*Alarms attribute*), 46
Alarm16 (*AlarmText attribute*), 44
Alarm17 (*Alarms attribute*), 46
Alarm17 (*AlarmText attribute*), 44, 67
Alarm18 (*Alarms attribute*), 46
Alarm18 (*AlarmText attribute*), 44
Alarm19 (*Alarms attribute*), 46
Alarm19 (*AlarmText attribute*), 44, 67
Alarm2 (*Alarms attribute*), 46
Alarm2 (*AlarmText attribute*), 44, 67
Alarm20 (*Alarms attribute*), 46
Alarm20 (*AlarmText attribute*), 44, 67
Alarm21 (*Alarms attribute*), 46
Alarm21 (*AlarmText attribute*), 44, 67
Alarm22 (*Alarms attribute*), 46
Alarm22 (*AlarmText attribute*), 44, 67
Alarm23 (*Alarms attribute*), 46
Alarm23 (*AlarmText attribute*), 44
Alarm24 (*Alarms attribute*), 46
Alarm24 (*AlarmText attribute*), 44
Alarm25 (*Alarms attribute*), 46
Alarm25 (*AlarmText attribute*), 44
Alarm26 (*Alarms attribute*), 46
Alarm26 (*AlarmText attribute*), 44
Alarm27 (*Alarms attribute*), 46
Alarm28 (*Alarms attribute*), 46
Alarm29 (*Alarms attribute*), 46
Alarm3 (*Alarms attribute*), 46
Alarm3 (*AlarmText attribute*), 44, 67
Alarm30 (*Alarms attribute*), 47
Alarm31 (*Alarms attribute*), 47
Alarm32 (*Alarms attribute*), 47
Alarm33 (*Alarms attribute*), 47
Alarm34 (*Alarms attribute*), 47
Alarm35 (*Alarms attribute*), 47
Alarm36 (*Alarms attribute*), 47
Alarm37 (*Alarms attribute*), 47
Alarm38 (*Alarms attribute*), 47
Alarm39 (*Alarms attribute*), 47
Alarm4 (*Alarms attribute*), 47
Alarm4 (*AlarmText attribute*), 44, 67
Alarm40 (*Alarms attribute*), 47
Alarm41 (*Alarms attribute*), 47
Alarm41 (*AlarmText attribute*), 44
Alarm42 (*Alarms attribute*), 47
Alarm42 (*AlarmText attribute*), 44
Alarm43 (*Alarms attribute*), 47
Alarm43 (*AlarmText attribute*), 44
Alarm44 (*Alarms attribute*), 47
Alarm44 (*AlarmText attribute*), 44
Alarm45 (*Alarms attribute*), 47
Alarm45 (*AlarmText attribute*), 44
Alarm46 (*Alarms attribute*), 47
Alarm46 (*AlarmText attribute*), 44
Alarm47 (*Alarms attribute*), 47
Alarm47 (*AlarmText attribute*), 44
Alarm48 (*Alarms attribute*), 47
Alarm48 (*AlarmText attribute*), 44
Alarm49 (*Alarms attribute*), 47
Alarm5 (*Alarms attribute*), 47
Alarm5 (*AlarmText attribute*), 44, 67
Alarm50 (*Alarms attribute*), 47
Alarm51 (*Alarms attribute*), 47
Alarm52 (*Alarms attribute*), 47
Alarm53 (*Alarms attribute*), 47
Alarm54 (*Alarms attribute*), 47
Alarm55 (*Alarms attribute*), 47
Alarm56 (*Alarms attribute*), 47
Alarm57 (*Alarms attribute*), 47
Alarm58 (*Alarms attribute*), 47
Alarm59 (*Alarms attribute*), 47
Alarm6 (*Alarms attribute*), 47
Alarm6 (*AlarmText attribute*), 44, 67
Alarm60 (*Alarms attribute*), 47
Alarm61 (*Alarms attribute*), 47
Alarm62 (*Alarms attribute*), 47
Alarm63 (*Alarms attribute*), 48
Alarm64 (*Alarms attribute*), 48
Alarm65 (*Alarms attribute*), 48
Alarm66 (*Alarms attribute*), 48
Alarm67 (*Alarms attribute*), 48
Alarm68 (*Alarms attribute*), 48
Alarm69 (*Alarms attribute*), 48
Alarm7 (*Alarms attribute*), 48
Alarm7 (*AlarmText attribute*), 44, 67
Alarm70 (*Alarms attribute*), 48

Alarm71 (*Alarms attribute*), 48
 Alarm72 (*Alarms attribute*), 48
 Alarm73 (*Alarms attribute*), 48
 Alarm74 (*Alarms attribute*), 48
 Alarm75 (*Alarms attribute*), 48
 Alarm76 (*Alarms attribute*), 48
 Alarm77 (*Alarms attribute*), 48
 Alarm78 (*Alarms attribute*), 48
 Alarm79 (*Alarms attribute*), 48
 Alarm8 (*Alarms attribute*), 48
 Alarm8 (*AlarmText attribute*), 44, 67
 Alarm80 (*Alarms attribute*), 48
 Alarm81 (*Alarms attribute*), 48
 Alarm82 (*Alarms attribute*), 48
 Alarm83 (*Alarms attribute*), 48
 Alarm84 (*Alarms attribute*), 48
 Alarm85 (*Alarms attribute*), 48
 Alarm86 (*Alarms attribute*), 48
 Alarm87 (*Alarms attribute*), 48
 Alarm88 (*Alarms attribute*), 48
 Alarm89 (*Alarms attribute*), 48
 Alarm9 (*Alarms attribute*), 48
 Alarm9 (*AlarmText attribute*), 44, 67
 Alarm90 (*Alarms attribute*), 48
 Alarm91 (*Alarms attribute*), 48
 Alarm92 (*Alarms attribute*), 48
 Alarm93 (*Alarms attribute*), 48
 Alarm94 (*Alarms attribute*), 48
 Alarm95 (*Alarms attribute*), 48
 Alarm96 (*Alarms attribute*), 49
 Alarm97 (*Alarms attribute*), 49
 Alarm98 (*Alarms attribute*), 49
 Alarm99 (*Alarms attribute*), 49
Alarms (*class in hvl_ccb.dev.supercube.constants*), 44
AlarmText (*class in hvl_ccb.dev.supercube.constants*), 43
AlarmText (*class in hvl_ccb.dev.supercube2015.constants*), 67
ALL (*TiePieOscilloscopeAutoResolutionModes attribute*), 92
all_measurements (*LuminoxMeasurementType attribute*), 180
all_measurements_types() (*LuminoxMeasurementType class method*), 180
amplitude (*TiePieGeneratorConfig property*), 88
ANALOG (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 138
analog_control (*FuGProbusVDIRegisters property*), 118
ANY (*LabJack.DeviceType attribute*), 130
ANY (*LJMCommunicationConfig.ConnectionType attribute*), 16
ANY (*LJMCommunicationConfig.DeviceType attribute*), 16
ANY (*TiePieOscilloscopeTriggerKind attribute*), 87
ARBITRARY (*TiePieGeneratorSignalType attribute*), 89
AsyncCommunicationProtocol (*class in hvl_ccb.comm.base*), 11
AsyncCommunicationProtocolConfig (*class in hvl_ccb.comm.base*), 13
ATM (*Pressure attribute*), 187
ATMOSPHERE (*Pressure attribute*), 187
attenuation (*CryLasAttenuator property*), 98
auto (*EarthingStickOperatingStatus attribute*), 51
AUTO (*LaudaProRp245eConfig.OperationModeEnum attribute*), 138
AUTO (*RTO1024.TriggerModes attribute*), 165
auto_install_mode (*T560 property*), 35
auto_install_mode (*T560Config attribute*), 36
auto_laser_on (*CryLasLaserConfig attribute*), 103
auto_resolution_mode (*TiePieOscilloscopeConfig property*), 92
AutoInstallMode (*class in hvl_ccb.dev.hIGHLAND_t560.base*), 32
AutoNumberNameEnum (*class in hvl_ccb.utils.enum*), 188

B

B (*SupercubeOpcEndpoint attribute*), 57, 74
BA (*NewportConfigCommands attribute*), 145
backlash_compensation (*NewportSMC100PPConfig attribute*), 152
backup_waveform() (*RTO1024 method*), 165
bar (*PfeifferTPG.PressureUnits attribute*), 158
BAR (*Pressure attribute*), 187
barometric_pressure (*LuminoxMeasurementType attribute*), 180
base_velocity (*NewportSMC100PPConfig attribute*), 152
BATH_TEMP (*LaudaProRp245eCommand attribute*), 136
baudrate (*CryLasAttenuatorSerialCommunicationConfig attribute*), 99
baudrate (*CryLasLaserSerialCommunicationConfig attribute*), 105
baudrate (*FuGSerialCommunicationConfig attribute*), 121
baudrate (*HeinzingerSerialCommunicationConfig attribute*), 127
baudrate (*LuminoxSerialCommunicationConfig attribute*), 181
baudrate (*MBW973SerialCommunicationConfig attribute*), 143
baudrate (*NewportSMC100PPSerialCommunicationConfig attribute*), 156
baudrate (*PfeifferTPGSerialCommunicationConfig attribute*), 163
baudrate (*SerialCommunicationConfig attribute*), 24
BH (*NewportConfigCommands attribute*), 145
board (*VisaCommunicationConfig attribute*), 31

BreakdownDetection (class *in* `hvl_ccb.dev.supercube.constants`), 49
BreakdownDetection (class *in* `hvl_ccb.dev.supercube2015.constants`), 67
bufsize (`TcpCommunicationConfig` attribute), 26
build_str() (`LaudaProRp245eCommand` method), 137
bytesize (`CryLasAttenuatorSerialCommunicationConfig` attribute), 99
bytesize (`CryLasLaserSerialCommunicationConfig` attribute), 106
bytesize (`FuGSerialCommunicationConfig` attribute), 121
bytesize (`HeinzingerSerialCommunicationConfig` attribute), 127
bytesize (`LuminoxSerialCommunicationConfig` attribute), 181
bytesize (`MBW973SerialCommunicationConfig` attribute), 143
bytesize (`NewportSMC100PPSerialCommunicationConfig` attribute), 156
bytesize (`PfeifferTPGSerialCommunicationConfig` attribute), 163
Bytesize (`SerialCommunicationConfig` attribute), 24
bytesize (`SerialCommunicationConfig` attribute), 24

C

C (`LabJack.TemperatureUnit` attribute), 130
C (`LabJack.ThermocoupleType` attribute), 130
C (`Temperature` attribute), 187
calibration_factor (`CryLasLaserConfig` attribute), 103
calibration_factor (`LEM4000S` attribute), 186
calibration_mode (`FuGProbusVDIRegisters` property), 118
cc_mode (`FuGProbusVDIRegisters` property), 118
CCBException, 192
cee16 (`GeneralSockets` attribute), 52, 70
CELSIUS (`Temperature` attribute), 187
ch_a (`T560` property), 35
ch_b (`T560` property), 35
ch_c (`T560` property), 35
ch_d (`T560` property), 35
channels_enabled (`TiePieOscilloscope` property), 91
check_for_error() (`NewportSMC100PPSerialCommunication` method), 154
check_generic_type() (*in module* `hvl_ccb.utils.typing`), 189
check_master_slave_config() (`PSI9000` method), 107
chunk_size (`VisaCommunicationConfig` attribute), 31
clean_amplitude() (`TiePieGeneratorConfig` method), 88
clean_auto_resolution_mode() (`TiePieOscilloscopeConfig` static method), 93
clean_coupling() (`TiePieOscilloscopeChannelConfig` static method), 85
clean_enabled() (`TiePieGeneratorConfig` static method), 88
clean_enabled() (`TiePieOscilloscopeChannelConfig` static method), 85
clean_frequency() (`TiePieGeneratorConfig` method), 88
clean_input_range() (`TiePieOscilloscopeChannelConfig` method), 85
clean_offset() (`TiePieGeneratorConfig` method), 88
clean_pre_sample_ratio() (`TiePieOscilloscopeConfig` method), 93
clean_probe_offset() (`TiePieOscilloscopeChannelConfig` method), 85
clean_record_length() (`TiePieOscilloscopeConfig` method), 93
clean_resolution() (`TiePieOscilloscopeConfig` static method), 93
clean_sample_frequency() (`TiePieOscilloscopeConfig` method), 93
clean_signal_type() (`TiePieGeneratorConfig` static method), 88
clean_trigger_enabled() (`TiePieOscilloscopeChannelConfig` static method), 85
clean_trigger_hysteresis() (`TiePieOscilloscopeChannelConfig` method), 85
clean_trigger_kind() (`TiePieOscilloscopeChannelConfig` static method), 85
clean_trigger_level() (`TiePieOscilloscopeChannelConfig` method), 85
clean_trigger_level_mode() (`TiePieOscilloscopeChannelConfig` static method), 85
clean_trigger_timeout() (`TiePieOscilloscopeConfig` method), 93
clean_values() (`AsyncCommunicationProtocolConfig` method), 13
clean_values() (`CryLasAttenuatorConfig` method), 98
clean_values() (`CryLasLaserConfig` method), 103
clean_values() (`EmptyConfig` method), 96, 191
clean_values() (`FuGConfig` method), 113
clean_values() (`HeinzingerConfig` method), 123
clean_values() (`ILS2TConfig` method), 175
clean_values() (`LaudaProRp245eConfig` method), 138
clean_values() (*in module* `LaudaProRp245eTcpCommunicationConfig` method), 140
clean_values() (`LJMCommunicationConfig` method), 17
clean_values() (`LuminoxConfig` method), 179
clean_values() (`MBW973Config` method), 142

clean_values() (*ModbusTcpCommunicationConfig method*), 19
clean_values() (*NewportSMC100PPConfig method*), 152
clean_values() (*OpcUaCommunicationConfig method*), 21
clean_values() (*PfeifferTPGConfig method*), 161
clean_values() (*PSI9000Config method*), 109
clean_values() (*SerialCommunicationConfig method*), 24
clean_values() (*SupercubeConfiguration method*), 41, 65
clean_values() (*T560Config method*), 36
clean_values() (*TcpCommunicationConfig method*), 26
clean_values() (*TechnixConfig method*), 81
clean_values() (*TelnetCommunicationConfig method*), 28
clean_values() (*TiePieDeviceConfig method*), 83
clean_values() (*VisaCommunicationConfig method*), 31
clean_waveform() (*TiePieGeneratorConfig method*), 88
close (*EarthingStickOperation attribute*), 51
close() (*CommunicationProtocol method*), 14
close() (*LaudaProRp245eTcpCommunication method*), 139
close() (*LJMCommunication method*), 15
close() (*ModbusTcpCommunication method*), 18
close() (*NullCommunicationProtocol method*), 14
close() (*OpcUaCommunication method*), 20
close() (*SerialCommunication method*), 23
close() (*Tcp method*), 26
close() (*TelnetCommunication method*), 28
close() (*VisaCommunication method*), 30
close_shutter() (*CryLasLaser method*), 101
CLOSED (*CryLasLaser.AnswersShutter attribute*), 100
CLOSED (*CryLasLaserShutterStatus attribute*), 106
closed (*DoorStatus attribute*), 49, 68
closed (*EarthingStickStatus attribute*), 52, 69
CMD_EXEC_ERROR (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_CC (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_CONFIGURATION (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_DISABLE (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
attribute), 154
CMD_NOT_ALLOWED_HOMING (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_MOVING (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_NOT_REFERENCED (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_PP (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMD_NOT_ALLOWED_READY (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
CMR (*PfeifferTPG.SensorTypes attribute*), 159
CODE_OR_ADDR_INVALID (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
collect_measurement_data() (*TiePieOscilloscope method*), 91
com (*SingleCommDevice property*), 97
COM_TIME_OUT (*LaudaProRp245eCommand attribute*), 136
com_time_out (*LaudaProRp245eConfig attribute*), 138
COM_TIMEOUT (*NewportSMC100PPSerialCommunication.ControllerErrors attribute*), 154
command (*LuminoxMeasurementType property*), 180
COMMAND (*TriggerMode attribute*), 34
command() (*FuGProbusIV method*), 116
communication_channel (*TechnixConfig attribute*), 81
CommunicationException, 14
CommunicationProtocol (*class in hvl_ccb.comm.base*), 14
config (*ConfigurationMixin property*), 190
CONFIG (*FuGProbusVRegisterGroups attribute*), 119
CONFIG (*NewportSMC100PP.StateMessages attribute*), 146
CONFIG (*NewportStates attribute*), 157
config_cls() (*AsyncCommunicationProtocol static config_cls() ConfigurationMixin static method*), 11
config_cls() (*ConfigurationMixin static method*), 190
config_cls() (*CryLasAttenuator static method*), 98
config_cls() (*CryLasAttenuatorSerialCommunication static method*), 99
config_cls() (*CryLasLaser static method*), 101
config_cls() (*CryLasLaserSerialCommunication static method*), 105
config_cls() (*Device static method*), 95
config_cls() (*FuGProbusIV static method*), 116
config_cls() (*FuGSerialCommunication static method*), 121
config_cls() (*HeinzingerDI static method*), 124

config_cls() (*HeinzingerSerialCommunication static method*), 126
config_cls() (*ILS2T static method*), 173
config_cls() (*ILS2TModbusTcpCommunication static method*), 176
config_cls() (*LaudaProRp245e static method*), 134
config_cls() (*LaudaProRp245eTcpCommunication static method*), 139
config_cls() (*LJMCommunication static method*), 15
config_cls() (*Luminox static method*), 178
config_cls() (*LuminoxSerialCommunication static method*), 181
config_cls() (*MBW973 static method*), 141
config_cls() (*MBW973SerialCommunication static method*), 143
config_cls() (*ModbusTcpCommunication static method*), 18
config_cls() (*NewportSMC100PP static method*), 146
config_cls() (*NewportSMC100PPSerialCommunication static method*), 154
config_cls() (*NullCommunicationProtocol static method*), 14
config_cls() (*OpcUaCommunication static method*), 20
config_cls() (*PfeifferTPG static method*), 159
config_cls() (*PfeifferTPGSerialCommunication static method*), 162
config_cls() (*PSI9000 static method*), 107
config_cls() (*PSI9000VisaCommunication static method*), 110
config_cls() (*RTO1024 static method*), 165
config_cls() (*RTO1024VisaCommunication static method*), 171
config_cls() (*SerialCommunication static method*), 23
config_cls() (*Supercube2015Base static method*), 62
config_cls() (*SupercubeAOpcUaCommunication static method*), 57, 75
config_cls() (*SupercubeBase static method*), 37
config_cls() (*SupercubeBOpcUaCommunication static method*), 60
config_cls() (*SupercubeOpcUaCommunication static method*), 42, 65
config_cls() (*SyncCommunicationProtocol static method*), 14
config_cls() (*T560 static method*), 35
config_cls() (*T560Communication static method*), 33
config_cls() (*Tcp static method*), 26
config_cls() (*Technix static method*), 80
config_cls() (*TechnixSerialCommunication static method*), 77
config_cls() (*TechnixTelnetCommunication static method*), 78
config_cls() (*TelnetCommunication static method*), 28
config_cls() (*TiePieOscilloscope static method*), 91
config_cls() (*VisaCommunication static method*), 30
config_cls() (*VisaDevice static method*), 184
config_gen() (*TiePieGeneratorMixin attribute*), 89
config_high_pulse() (*LabJack method*), 131
config_i2c() (*TiePieI2CHostMixin attribute*), 90
config_osc() (*TiePieOscilloscope attribute*), 91
config_osc_channel_dict (*TiePieOscilloscope attribute*), 91
config_status() (*FuG property*), 111
configdataclass() (*in module hvl_ccb.configuration*), 191
configuration_save_json() (*ConfigurationMixin method*), 190
ConfigurationMixin (*class in hvl_ccb.configuration*), 190
ConfigurationValueWarning, 190
connection_type (*LJMCommunicationConfig attribute*), 17
CONT_MODE (*LaudaProRp245eCommand attribute*), 137
contact_range() (*GeneralSupport class method*), 52
continue_ramp() (*LaudaProRp245e method*), 135
control_mode (*LaudaProRp245eConfig attribute*), 138
CONVERSION (*LEM4000S attribute*), 185
convert() (*LEM4000S method*), 186
convert() (*LMT70A method*), 186
convert() (*Sensor method*), 186
convert() (*Unit class method*), 187
ConvertibleTypes (*in module hvl_ccb.utils.typing*), 189
COOLOFF (*LaudaProRp245eConfig.OperationModeEnum attribute*), 138
COOLON (*LaudaProRp245eConfig.OperationModeEnum attribute*), 138
coupling (*TiePieOscilloscopeChannelConfig property*), 85
CR (*FuGTerminators attribute*), 122
create_serial_port() (*SerialCommunicationConfig method*), 24
create_telnet() (*TelnetCommunicationConfig method*), 28
CRLF (*FuGTerminators attribute*), 122
CryLasAttenuator (*class in hvl_ccb.dev.crylas*), 98
CryLasAttenuatorConfig (*class in hvl_ccb.dev.crylas*), 98
CryLasAttenuatorError, 99
CryLasAttenuatorSerialCommunication (*class in hvl_ccb.dev.crylas*), 99
CryLasAttenuatorSerialCommunicationConfig (*class in hvl_ccb.dev.crylas*), 99
CryLasLaser (*class in hvl_ccb.dev.crylas*), 100
CryLasLaser.AnswersShutter (*class in hvl_ccb.dev.crylas*), 100
CryLasLaser.AnswersStatus (*class in hvl_ccb.dev.crylas*), 100

CryLasLaser.LaserStatus	(class in hvl_ccb.dev.crylas), 101	default_com_cls()	(NewportSMC100PP static method), 146
CryLasLaser.RepetitionRates	(class in hvl_ccb.dev.crylas), 101	default_com_cls()	(PfeifferTPG static method), 159
CryLasLaserConfig (class in hvl_ccb.dev.crylas)	, 103	default_com_cls()	(PSI9000 static method), 107
CryLasLaserError	, 104	default_com_cls()	(RTO1024 static method), 165
CryLasLaserNotReadyError	, 104	default_com_cls()	(SingleCommDevice static method), 97
CryLasLaserPoller (class in hvl_ccb.dev.crylas)	, 104	default_com_cls()	(Supercube2015Base static method), 62
CryLasLaserSerialCommunication	(class in hvl_ccb.dev.crylas), 104	default_com_cls()	(Supercube2015WithFU static method), 74
CryLasLaserSerialCommunicationConfig (class in hvl_ccb.dev.crylas), 105		default_com_cls()	(SupercubeB static method), 60
CryLasLaserShutterStatus	(class in hvl_ccb.dev.crylas), 106	default_com_cls()	(SupercubeBase static method), 37
current (FuG property)	, 111	default_com_cls()	(SupercubeWithFU static method), 58
CURRENT (FuGProbusIVCommands attribute)	, 116	default_com_cls()	(T560 static method), 35
CURRENT (FuGReadbackChannels attribute)	, 120	default_com_cls()	(Technix method), 80
current (Technix property)	, 80	default_com_cls()	(TiePieOscilloscope static method), 92
current_lower_limit (PSI9000Config attribute)	, 109	default_com_cls()	(VisaDevice static method), 184
current_monitor (FuG property)	, 112	DEFAULT_IO_SCANNING_CONTROL_VALUES (ILS2T attribute)	, 172
current_primary (Power attribute)	, 55, 72	default_n_attempts_read_text_nonempty (AsyncCommunicationProtocolConfig attribute)	, 13
current_upper_limit (PSI9000Config attribute)	, 109	default_n_attempts_read_text_nonempty (FuGSerialCommunicationConfig attribute)	, 121
CurrentPosition	(NewportSMC100PPConfig.HomeSearch attribute), 152	default_n_attempts_read_text_nonempty (HeinzingerSerialCommunicationConfig attribute)	, 127
cv_mode (FuGProbusVDIRegisters property)	, 118	default_number_of_recordings (HeinzingerConfig attribute)	, 123
D		Device (class in hvl_ccb.dev.base)	, 95
datachange_notification() (OpcUaSubHandler method)	, 22	DEVICE_TYPE (LaudaProRp245eCommand attribute)	, 137
datachange_notification() (SupercubeSubscriptionHandler method)	, 43, 66	device_type (LJMCommunicationConfig attribute)	, 17
date_of_manufacture (LuminoxMeasurementType attribute)	, 180	DeviceException	, 95
DC (TiePieGeneratorSignalType attribute)	, 89	DeviceExistingException	, 95
DC_DoubleStage_280kV (PowerSetup attribute)	, 55, 72	DeviceFailuresException	, 95
DC_SingleStage_140kV (PowerSetup attribute)	, 56, 73	devices_failed_start (DeviceSequenceMixin attribute)	, 96
DC_VOLTAGE_TOO_LOW	(NewportSMC100PP.MotorErrors attribute), 145	devices_failed_start (ExperimentManager attribute)	, 193
DCA (TiePieOscilloscopeChannelCoupling attribute)	, 86	devices_failed_stop (DeviceSequenceMixin attribute)	, 96
DCV (TiePieOscilloscopeChannelCoupling attribute)	, 86	devices_failed_stop (ExperimentManager attribute)	, 193
default_com_cls() (CryLasAttenuator static method)	, 98	DeviceSequenceMixin (class in hvl_ccb.dev.base)	, 95
default_com_cls() (CryLasLaser static method)	, 101	di (FuG property)	, 112
default_com_cls() (FuGProbusIV static method)	, 116	digital_control (FuGProbusVDIRegisters property)	, 118
default_com_cls() (HeinzingerDI static method)	, 124	DIOChannel (LabJack attribute)	, 129
default_com_cls() (ILS2T static method)	, 173	DISABLE (NewportStates attribute)	, 157
default_com_cls() (LabJack static method)	, 131	disable() (ILS2T method)	, 173
default_com_cls() (LaudaProRp245e static method)	, 135	DISABLE_FROM_JOGGING	(New-
default_com_cls() (Luminox static method)	, 179		
default_com_cls() (MBW973 static method)	, 141		

<i>portSMC100PP.StateMessages</i>	<i>attribute), E505 (FuGErrorcodes attribute), 115</i>
146	<i>E6 (FuGErrorcodes attribute), 115</i>
DISABLE_FROM_MOVING	<i>(New-attribute), E666 (FuGErrorcodes attribute), 115</i>
<i>portSMC100PP.StateMessages</i>	<i>E7 (FuGErrorcodes attribute), 115</i>
146	<i>E8 (FuGErrorcodes attribute), 115</i>
DISABLE_FROM_READY	<i>(New-attribute), E9 (FuGErrorcodes attribute), 115</i>
<i>portSMC100PP.StateMessages</i>	<i>EarthingRod (class in hvl_ccb.dev.supercube.constants), 49</i>
146	<i>EarthingRodStatus (class in hvl_ccb.dev.supercube.constants), 50</i>
disable_pulses() (LabJack method), 131	<i>EarthingStick (class in hvl_ccb.dev.supercube.constants), 50</i>
DISABLED (TiePieOscilloscopeAutoResolutionModes attribute), 92	<i>EarthingStick (class in hvl_ccb.dev.supercube2015.constants), 68</i>
DisableEspStageCheck	<i>EarthingStickMeta (class in hvl_ccb.dev.supercube.constants), 51</i>
<i>portSMC100PPConfig.EspStageConfig attribute), 152</i>	<i>EarthingStickOperatingStatus (class in hvl_ccb.dev.supercube.constants), 51</i>
disarm_trigger() (T560 method), 36	<i>EarthingStickOperation (class in hvl_ccb.dev.supercube.constants), 51</i>
DISPLACEMENT_OUT_OF_LIMIT	<i>EarthingStickStatus (class in hvl_ccb.dev.supercube.constants), 51</i>
<i>portSMC100PPSerialCommunication.Controller attribute), 154</i>	<i>EarthingStickStatus (class in hvl_ccb.dev.supercube2015.constants), 69</i>
display_message_board() (SupercubeBase method), 38	EEPROM_ACCESS_ERROR (New-portSMC100PPSerialCommunication.ControllerErrors attribute), 154
display_status_board() (SupercubeBase method), 38	EIGHT (HeinzingerConfig.RecordingsEnum attribute), 123
do_iostressing_write() (ILS2T method), 173	EIGHT_BIT (TiePieOscilloscopeResolution attribute), 93
Door (class in hvl_ccb.dev.supercube.constants), 49	EIGHT_HUNDRED_MILLI_VOLT (TiePieOscilloscopeRange attribute), 86
DoorStatus (class in hvl_ccb.dev.supercube.constants), 49	EIGHT_VOLT (TiePieOscilloscopeRange attribute), 86
DoorStatus (class in hvl_ccb.dev.supercube2015.constants), 68	EIGHTBITS (SerialCommunicationBytesize attribute), 23
E	EIGHTY_VOLT (TiePieOscilloscopeRange attribute), 86
E (LabJack.ThermocoupleType attribute), 130	EmptyConfig (class in hvl_ccb.configuration), 191
E0 (FuGErrorcodes attribute), 114	EmptyConfig (class in hvl_ccb.dev.base), 96
E1 (FuGErrorcodes attribute), 114	enable() (ILS2T method), 173
E10 (FuGErrorcodes attribute), 114	enable_clock() (LabJack method), 131
E100 (FuGErrorcodes attribute), 114	enabled (TiePieGeneratorConfig property), 88
E106 (FuGErrorcodes attribute), 114	enabled (TiePieOscilloscopeChannelConfig property), 85
E11 (FuGErrorcodes attribute), 114	EnableEspStageCheck (New-portSMC100PPConfig.EspStageConfig attribute), 152
E115 (FuGErrorcodes attribute), 114	encoding (AsyncCommunicationProtocolConfig attribute), 13
E12 (FuGErrorcodes attribute), 114	encoding (NewportSMC100PPSerialCommunicationConfig attribute), 156
E125 (FuGErrorcodes attribute), 114	encoding_error_handling (AsyncCommunicationProtocolConfig attribute), 13
E13 (FuGErrorcodes attribute), 114	encoding_error_handling (New-portSMC100PPSerialCommunicationConfig attribute), 156
E135 (FuGErrorcodes attribute), 114	
E14 (FuGErrorcodes attribute), 114	
E145 (FuGErrorcodes attribute), 114	
E15 (FuGErrorcodes attribute), 114	
E155 (FuGErrorcodes attribute), 114	
E16 (FuGErrorcodes attribute), 115	
E165 (FuGErrorcodes attribute), 115	
E2 (FuGErrorcodes attribute), 115	
E206 (FuGErrorcodes attribute), 115	
E306 (FuGErrorcodes attribute), 115	
E4 (FuGErrorcodes attribute), 115	
E5 (FuGErrorcodes attribute), 115	
E504 (FuGErrorcodes attribute), 115	

F

F (*LabJack.TemperatureUnit attribute*), 130
 F (*Temperature attribute*), 187
FAHRENHEIT (*Temperature attribute*), 187
 failures (*DeviceFailuresException attribute*), 95
FALLING (*TiePieOscilloscopeTriggerKind attribute*), 87
FAST (*ILS2T.Ref16Jog attribute*), 172
 file_copy() (*RTO1024 method*), 165
 finish() (*ExperimentManager method*), 193
FINISHED (*ExperimentStatus attribute*), 193
FINISHING (*ExperimentStatus attribute*), 193
 fire_trigger() (*T560 method*), 36
FIRMWARE (*FuGReadbackChannels attribute*), 120
FIVE_MHZ (*LabJack.ClockFrequency attribute*), 129
FIVEBITS (*SerialCommunicationBytesize attribute*), 23
FLT_INFO (*ILS2TRegAddr attribute*), 177
FLT_MEM_DEL (*ILS2TRegAddr attribute*), 177
FLT_MEM_RESET (*ILS2TRegAddr attribute*), 177
FOLLOWING_ERROR (*NewportSMC100PP.MotorErrors attribute*), 145
FOLLOWRAMP (*FuGRampModes attribute*), 120
force_trigger() (*TiePieOscilloscope method*), 92
 force_value() (*AsyncCommunicationProtocolConfig method*), 13
 force_value() (*CryLasAttenuatorConfig method*), 98
 force_value() (*CryLasAttenuatorSerialCommunicationConfig method*), 99
 force_value() (*CryLasLaserConfig method*), 104
 force_value() (*CryLasLaserSerialCommunicationConfig method*), 106
 force_value() (*EmptyConfig method*), 96, 191
 force_value() (*FuGConfig method*), 113
 force_value() (*FuGSerialCommunicationConfig method*), 121
 force_value() (*HeinzingerConfig method*), 123
 force_value() (*HeinzingerSerialCommunicationConfig method*), 127
 force_value() (*ILS2TConfig method*), 175
 force_value() (*ILS2TModbusTcpCommunicationConfig method*), 176
 force_value() (*LaudaProRp245eConfig method*), 138
 force_value() (*LaudaProRp245eTcpCommunicationConfig method*), 140
 force_value() (*LJMCommunicationConfig method*), 17
 force_value() (*LuminoxConfig method*), 179
 force_value() (*LuminoxSerialCommunicationConfig method*), 182
 force_value() (*MBW973Config method*), 142
 force_value() (*MBW973SerialCommunicationConfig method*), 143
 force_value() (*ModbusTcpCommunicationConfig method*), 19

force_value() (*NewportSMC100PPConfig method*), 152
force_value() (*NewportSMC100PPSerialCommunicationConfig method*), 156
force_value() (*OpcUaCommunicationConfig method*), 21
force_value() (*PfeifferTPGConfig method*), 161
force_value() (*PfeifferTPGSerialCommunicationConfig method*), 163
force_value() (*PSI9000Config method*), 109
force_value() (*PSI9000VisaCommunicationConfig method*), 110
force_value() (*RTO1024Config method*), 170
force_value() (*RTO1024VisaCommunicationConfig method*), 171
force_value() (*SerialCommunicationConfig method*), 24
force_value() (*SupercubeAOpcUaConfiguration method*), 58, 76
force_value() (*SupercubeBOpcUaConfiguration method*), 60
force_value() (*SupercubeConfiguration method*), 41, 65
force_value() (*SupercubeOpcUaCommunicationConfig method*), 42, 65
force_value() (*T560CommunicationConfig method*), 33
force_value() (*T560Config method*), 36
force_value() (*TcpCommunicationConfig method*), 26
force_value() (*TechnixConfig method*), 81
force_value() (*TechnixSerialCommunicationConfig method*), 78
force_value() (*TechnixTelnetCommunicationConfig method*), 79
force_value() (*TelnetCommunicationConfig method*), 28
force_value() (*TiePieDeviceConfig method*), 83
force_value() (*VisaCommunicationConfig method*), 31
force_value() (*VisaDeviceConfig method*), 184
FORTY_MHZ (*LabJack.ClockFrequency attribute*), 129
FORTY_VOLT (*TiePieOscilloscopeRange attribute*), 86
FOUR (*HeinzingerConfig.RecordingsEnum attribute*), 123
FOUR_HUNDRED_MILLI_VOLT (*TiePieOscilloscopeRange attribute*), 86
FOUR_VOLT (*TiePieOscilloscopeRange attribute*), 86
FOURTEEN_BIT (*TiePieOscilloscopeResolution attribute*), 93
FREERUN (*RTO1024.TriggerModes attribute*), 165
frequency (*Power attribute*), 55, 72
frequency (*T560 property*), 36
frequency (*TiePieGeneratorConfig property*), 88
FRM (*NewportConfigCommands attribute*), 145
from_json() (*ConfigurationMixin class method*), 190
FRS (*NewportConfigCommands attribute*), 145
fso_reset() (*Supercube2015WithFU method*), 74
fso_reset() (*SupercubeWithFU method*), 58
FuG (*class in hvl_ccb.dev.fug*), 111
FuGConfig (*class in hvl_ccb.dev.fug*), 113
FuGDigitalVal (*class in hvl_ccb.dev.fug*), 113
FuGError, 114
FuGErrorcodes (*class in hvl_ccb.dev.fug*), 114
FuGMonitorModes (*class in hvl_ccb.dev.fug*), 115
FuGPolarities (*class in hvl_ccb.dev.fug*), 116
FuGProbusIV (*class in hvl_ccb.dev.fug*), 116
FuGProbusIVCommands (*class in hvl_ccb.dev.fug*), 116
FuGProbusV (*class in hvl_ccb.dev.fug*), 117
FuGProbusVConfigRegisters (*class in hvl_ccb.dev.fug*), 117
FuGProbusVDIRegisters (*class in hvl_ccb.dev.fug*), 118
FuGProbusVDORegisters (*class in hvl_ccb.dev.fug*), 118
FuGProbusVMonitorRegisters (*class in hvl_ccb.dev.fug*), 119
FuGProbusVRegisterGroups (*class in hvl_ccb.dev.fug*), 119
FuGProbusVSetRegisters (*class in hvl_ccb.dev.fug*), 119
FuGRampModes (*class in hvl_ccb.dev.fug*), 120
FuGReadbackChannels (*class in hvl_ccb.dev.fug*), 120
FuGSerialCommunication (*class in hvl_ccb.dev.fug*), 120
FuGSerialCommunicationConfig (*class in hvl_ccb.dev.fug*), 121
FuGTerminators (*class in hvl_ccb.dev.fug*), 122

G

gate_mode (*T560 property*), 36
gate_polarity (*T560 property*), 36
GateMode (*class in hvl_ccb.dev.highland_t560.base*), 33
GeneralSockets (*class in hvl_ccb.dev.supercube.constants*), 52
GeneralSockets (*class in hvl_ccb.dev.supercube2015.constants*), 70
GeneralSupport (*class in hvl_ccb.dev.supercube.constants*), 52
GeneralSupport (*class in hvl_ccb.dev.supercube2015.constants*), 70
GeneralSupportMeta (*class in hvl_ccb.dev.supercube.constants*), 54
GENERATOR (*TiePieDeviceType attribute*), 84
generator_start() (*TiePieGeneratorMixin method*), 89
generator_stop() (*TiePieGeneratorMixin method*), 89
get() (*AlarmText class method*), 44, 67
get() (*MeasurementsDividerRatio class method*), 71

get() (*MeasurementsScaledInput* class method), 72
 get_acceleration() (*NewportSMC100PP* method), 147
 get_acquire_length() (*RTO1024* method), 165
 get_ain() (*LabJack* method), 131
 get_bath_temp() (*LaudaProRp245e* method), 135
 get_by_p_id() (*LabJack.DeviceType* class method), 130
 get_by_p_id() (*LJMCommunicationConfig.DeviceType* class method), 16
 get_cal_current_source() (*LabJack* method), 131
 get_cee16_socket() (*Supercube2015Base* method), 62
 get_cee16_socket() (*SupercubeBase* method), 38
 get_channel_offset() (*RTO1024* method), 165
 get_channel_position() (*RTO1024* method), 166
 get_channel_range() (*RTO1024* method), 166
 get_channel_scale() (*RTO1024* method), 166
 get_channel_state() (*RTO1024* method), 166
 get_clock() (*LabJack* method), 131
 get_controller_information() (*NewportSMC100PP* method), 147
 get_current() (*HeinzingerDI* method), 124
 get_dc_volt() (*ILS2T* method), 174
 get_device() (*DeviceSequenceMixin* method), 96
 get_device_by_serial_number() (in *module hvl_ccb.dev.tiepie.base*), 84
 get_device_type() (*LaudaProRp245e* method), 135
 get_devices() (*DeviceSequenceMixin* method), 96
 get_digital_input() (*LabJack* method), 131
 get_door_status() (*Supercube2015Base* method), 62
 get_door_status() (*SupercubeBase* method), 38
 get_earthing_manual() (*Supercube2015Base* method), 62
 get_earthing_rod_status() (*SupercubeBase* method), 38
 get_earthing_status() (*Supercube2015Base* method), 62
 get_earthing_stick_manual() (*SupercubeBase* method), 38
 get_earthing_stick_operating_status() (*SupercubeBase* method), 38
 get_earthing_stick_status() (*SupercubeBase* method), 38
 get_error_code() (*ILS2T* method), 174
 get_error_queue() (*VisaDevice* method), 184
 get_frequency() (*Supercube2015WithFU* method), 75
 get_frequency() (*SupercubeWithFU* method), 58
 get_fso_active() (*Supercube2015WithFU* method), 75
 get_fso_active() (*SupercubeWithFU* method), 58
 get_full_scale_mbar() (*PfeifferTPG* method), 159
 get_full_scale_unitless() (*PfeifferTPG* method), 159

get_identification() (*VisaDevice* method), 184
 get_interface_version() (*HeinzingerDI* method), 124
 get_max_voltage() (*Supercube2015WithFU* method), 75
 get_max_voltage() (*SupercubeWithFU* method), 58
 get_measurement_ratio() (*Supercube2015Base* method), 62
 get_measurement_ratio() (*SupercubeBase* method), 38
 get_measurement_voltage() (*Supercube2015Base* method), 62
 get_measurement_voltage() (*SupercubeBase* method), 38
 get_motor_configuration() (*NewportSMC100PP* method), 147
 get_move_duration() (*NewportSMC100PP* method), 147
 get_negative_software_limit() (*NewportSMC100PP* method), 148
 get_number_of_recordings() (*HeinzingerDI* method), 124
 get_output() (*PSI9000* method), 107
 get_position() (*ILS2T* method), 174
 get_position() (*NewportSMC100PP* method), 148
 get_positive_software_limit() (*NewportSMC100PP* method), 148
 get_power_setup() (*Supercube2015WithFU* method), 75
 get_power_setup() (*SupercubeWithFU* method), 58
 get_primary_current() (*Supercube2015WithFU* method), 75
 get_primary_current() (*SupercubeWithFU* method), 59
 get_primary_voltage() (*Supercube2015WithFU* method), 75
 get_primary_voltage() (*SupercubeWithFU* method), 59
 get_product_id() (*LabJack* method), 131
 get_product_name() (*LabJack* method), 131
 get_product_type() (*LabJack* method), 132
 get_pulse_energy_and_rate() (*CryLasLaser* method), 101
 get_reference_point() (*RTO1024* method), 166
 get_register() (*FuGProbusV* method), 117
 get_repetitions() (*RTO1024* method), 166
 get_sbus_rh() (*LabJack* method), 132
 get_sbus_temp() (*LabJack* method), 132
 get_serial_number() (*HeinzingerDI* method), 124
 get_serial_number() (*LabJack* method), 132
 get_state() (*NewportSMC100PP* method), 148
 get_status() (*ILS2T* method), 174
 get_status() (*Supercube2015Base* method), 63
 get_status() (*SupercubeBase* method), 39

```

get_support_input() (Supercube2015Base method), 63
get_support_input() (SupercubeBase method), 39
get_support_output() (Supercube2015Base method), 63
get_support_output() (SupercubeBase method), 39
get_system_lock() (PSI9000 method), 107
get_t13_socket() (Supercube2015Base method), 63
get_t13_socket() (SupercubeBase method), 39
get_target_voltage() (Supercube2015WithFU method), 75
get_target_voltage() (SupercubeWithFU method), 59
get_temperature() (ILS2T method), 174
get_timestamps() (RTO1024 method), 166
get_ui_lower_limits() (PSI9000 method), 108
get_uip_upper_limits() (PSI9000 method), 108
get_voltage() (HeinzingerDI method), 124
get_voltage_current_setpoint() (PSI9000 method), 108
go_home() (NewportSMC100PP method), 149
go_to_configuration() (NewportSMC100PP method), 149
GreenNotReady (SafetyStatus attribute), 56, 73
GreenReady (SafetyStatus attribute), 56, 73

H
HARDWARE (CryLasLaser.RepetitionRates attribute), 101
has_safe_ground (TiePieOscilloscopeChannelConfig property), 85
HEAD (CryLasLaser.AnswersStatus attribute), 100
HeinzingerConfig (class in hvl_ccb.dev.heinzinger), 122
HeinzingerConfig.RecordingsEnum (class in hvl_ccb.dev.heinzinger), 123
HeinzingerDI (class in hvl_ccb.dev.heinzinger), 123
HeinzingerDI.OutputStatus (class in hvl_ccb.dev.heinzinger), 123
HeinzingerPNC (class in hvl_ccb.dev.heinzinger), 125
HeinzingerPNC.UnitCurrent (class in hvl_ccb.dev.heinzinger), 125
HeinzingerPNC.UnitVoltage (class in hvl_ccb.dev.heinzinger), 125
HeinzingerPNCDeviceNotRecognizedException, 126
HeinzingerPNCError, 126
HeinzingerPNCMaxCurrentExceededException, 126
HeinzingerPNCMaxVoltageExceededException, 126
HeinzingerSerialCommunication (class in hvl_ccb.dev.heinzinger), 126
HeinzingerSerialCommunicationConfig (class in hvl_ccb.dev.heinzinger), 126
HIGH (LabJack.DIOStatus attribute), 130
high_resolution (FuGProbusVSetRegisters property), 119
home_search_polling_interval (NewportSMC100PPConfig attribute), 153
home_search_timeout (NewportSMC100PPConfig attribute), 153
home_search_type (NewportSMC100PPConfig attribute), 153
home_search_velocity (NewportSMC100PPConfig attribute), 153
HOME_STARTED (NewportSMC100PPSerialCommunication.ControllerError attribute), 154
HomeSwitch (NewportSMC100PPConfig.HomeSearch attribute), 152
HomeSwitch_and_Index (NewportSMC100PPConfig.HomeSearch attribute), 152
HOMING (NewportStates attribute), 157
HOMING_FROM_RS232 (NewportSMC100PP.StateMessages attribute), 146
HOMING_FROM_SMC (NewportSMC100PP.StateMessages attribute), 146
HOMING_TIMEOUT (NewportSMC100PP.MotorErrors attribute), 145
horn (Safety attribute), 73
horn() (Supercube2015Base method), 63
host (ModbusTcpCommunicationConfig attribute), 19
host (OpcUaCommunicationConfig attribute), 21
host (TcpCommunicationConfig attribute), 27
host (TelnetCommunicationConfig attribute), 28
host (VisaCommunicationConfig attribute), 31
hPascal (PfeifferTPG.PressureUnits attribute), 158
HT (NewportConfigCommands attribute), 145
hvl_ccb
    module, 194
hvl_ccb.comm
    module, 32
hvl_ccb.comm.base
    module, 11
hvl_ccb.comm.labjack_ljm
    module, 15
hvl_ccb.comm.modbus_tcp
    module, 18
hvl_ccb.comm.opc
    module, 20
hvl_ccb.comm.serial
    module, 22
hvl_ccb.comm.tcp
    module, 26
hvl_ccb.comm.telnet
    module, 27
hvl_ccb.comm.visa
    module, 29

```

hvl_ccb.configuration
 module, 190
 hvl_ccb.dev
 module, 185
 hvl_ccb.dev.base
 module, 95
 hvl_ccb.dev.crylas
 module, 97
 hvl_ccb.dev.ea_psi9000
 module, 107
 hvl_ccb.dev.fug
 module, 111
 hvl_ccb.dev.heinzinger
 module, 122
 hvl_ccb.dev.highland_t560
 module, 37
 hvl_ccb.dev.highland_t560.base
 module, 32
 hvl_ccb.dev.highland_t560.channel
 module, 34
 hvl_ccb.dev.highland_t560.device
 module, 35
 hvl_ccb.dev.labjack
 module, 128
 hvl_ccb.dev.lauda
 module, 134
 hvl_ccb.dev.mbw973
 module, 141
 hvl_ccb.dev.newport
 module, 144
 hvl_ccb.dev.pfeiffer_tpg
 module, 158
 hvl_ccb.dev.rs_rto1024
 module, 164
 hvl_ccb.dev.se_ilst2t
 module, 172
 hvl_ccb.dev.sst_luminox
 module, 178
 hvl_ccb.dev.supercube
 module, 61
 hvl_ccb.dev.supercube.base
 module, 37
 hvl_ccb.dev.supercube.constants
 module, 43
 hvl_ccb.dev.supercube.typ_a
 module, 57
 hvl_ccb.dev.supercube.typ_b
 module, 60
 hvl_ccb.dev.supercube2015
 module, 76
 hvl_ccb.dev.supercube2015.base
 module, 61
 hvl_ccb.dev.supercube2015.constants
 module, 67

hvl_ccb.dev.supercube2015.typ_a
 module, 74
 hvl_ccb.dev.technix
 module, 82
 hvl_ccb.dev.technix.base
 module, 77
 hvl_ccb.dev.technix.device
 module, 80
 hvl_ccb.dev.tiepie
 module, 94
 hvl_ccb.dev.tiepie.base
 module, 83
 hvl_ccb.dev.tiepie.channel
 module, 85
 hvl_ccb.dev.tiepie.device
 module, 87
 hvl_ccb.dev.tiepie.generator
 module, 88
 hvl_ccb.dev.tiepie.i2c
 module, 90
 hvl_ccb.dev.tiepie.oscilloscope
 module, 91
 hvl_ccb.dev.tiepie.utils
 module, 93
 hvl_ccb.dev.utils
 module, 182
 hvl_ccb.dev.visa
 module, 183
 hvl_ccb.exception
 module, 192
 hvl_ccb.experiment_manager
 module, 192
 hvl_ccb.utils
 module, 190
 hvl_ccb.utils.conversion_sensor
 module, 185
 hvl_ccb.utils.conversion_unit
 module, 186
 hvl_ccb.utils.enum
 module, 188
 hvl_ccb.utils.typing
 module, 189
 hvl_ccb.utils.validation
 module, 189

hysteresis_compensation *(New-portSMC100PPConfig attribute)*, 153

|

I2C (*TiePieDeviceType attribute*), 84
 ID (*FuGProbusIVCommands attribute*), 117
 Identification_error (*PfeifferTPG.SensorStatus attribute*), 158
 identifier (*LJMCommunicationConfig attribute*), 17
 identify_device() (*FuG method*), 112

identify_device() (*HeinzingerPNC method*), 126
identify_sensors() (*PfeifferTPG method*), 159
IKR (*PfeifferTPG.SensorTypes attribute*), 159
IKR11 (*PfeifferTPG.SensorTypes attribute*), 159
IKR9 (*PfeifferTPG.SensorTypes attribute*), 159
ILS2T (*class in hvl_ccb.dev.se_ilst2t*), 172
ILS2T.ActionsPtp (*class in hvl_ccb.dev.se_ilst2t*), 172
ILS2T.Mode (*class in hvl_ccb.dev.se_ilst2t*), 172
ILS2T.Ref16Jog (*class in hvl_ccb.dev.se_ilst2t*), 172
ILS2T.State (*class in hvl_ccb.dev.se_ilst2t*), 173
ILS2TConfig (*class in hvl_ccb.dev.se_ilst2t*), 175
ILS2TException, 176
ILS2TModbusTcpCommunication (*class in hvl_ccb.dev.se_ilst2t*), 176
ILS2TModbusTcpCommunicationConfig (*class in hvl_ccb.dev.se_ilst2t*), 176
ILS2TRegAddr (*class in hvl_ccb.dev.se_ilst2t*), 177
ILS2TRegDatatype (*class in hvl_ccb.dev.se_ilst2t*), 177
IMMEDIATELY (*FuGRampModes attribute*), 120
IMR (*PfeifferTPG.SensorTypes attribute*), 159
in_1_1 (*GeneralSupport attribute*), 52, 70
in_1_2 (*GeneralSupport attribute*), 52, 70
in_2_1 (*GeneralSupport attribute*), 53, 70
in_2_2 (*GeneralSupport attribute*), 53, 70
in_3_1 (*GeneralSupport attribute*), 53, 70
in_3_2 (*GeneralSupport attribute*), 53, 70
in_4_1 (*GeneralSupport attribute*), 53, 70
in_4_2 (*GeneralSupport attribute*), 53, 70
in_5_1 (*GeneralSupport attribute*), 53, 70
in_5_2 (*GeneralSupport attribute*), 53, 70
in_6_1 (*GeneralSupport attribute*), 53, 70
in_6_2 (*GeneralSupport attribute*), 53, 70
INACTIVE (*CryLasLaser.AnswersStatus attribute*), 100
inactive (*DoorStatus attribute*), 49, 68
inactive (*EarthingStickStatus attribute*), 52, 69
inhibit (*Technix property*), 80
init_attenuation (*CryLasAttenuatorConfig attribute*), 98
init_monitored_nodes() (*OpcUaCommunication method*), 20
init_shutter_status (*CryLasLaserConfig attribute*), 104
initialize() (*NewportSMC100PP method*), 149
INITIALIZED (*ExperimentStatus attribute*), 193
INITIALIZING (*ExperimentStatus attribute*), 193
Initializing (*SafetyStatus attribute*), 56, 73
INPUT (*FuGProbusVRegisterGroups attribute*), 119
INPUT (*GateMode attribute*), 33
input() (*GeneralSupport class method*), 53, 70
input_1 (*MeasurementsDividerRatio attribute*), 54, 71
input_1 (*MeasurementsScaledInput attribute*), 54, 72
input_2 (*MeasurementsDividerRatio attribute*), 54
input_2 (*MeasurementsScaledInput attribute*), 54, 72
input_3 (*MeasurementsDividerRatio attribute*), 54
input_3 (*MeasurementsScaledInput attribute*), 54, 72
input_4 (*MeasurementsDividerRatio attribute*), 54
input_4 (*MeasurementsScaledInput attribute*), 54, 72
input_range (*TiePieOscilloscopeChannelConfig property*), 86
INSTALL (*AutoInstallMode attribute*), 33
INT32 (*ILS2TRegDatatype attribute*), 178
INT_SYNTHESIZER (*TriggerMode attribute*), 34
interface_type (*PSI9000VisaCommunicationConfig attribute*), 110
interface_type (*RTO1024VisaCommunicationConfig attribute*), 171
interface_type (*VisaCommunicationConfig attribute*), 31
internal (*LabJack.CjcType attribute*), 129
INTERNAL (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 138
Internal (*PowerSetup attribute*), 56, 73
InvalidSupercubeStatusError, 61
IO_SCANNING (*ILS2TRegAddr attribute*), 177
IoScanningModeValueError, 178
is_configdataclass (*AsyncCommunicationProtocolConfig attribute*), 13
is_configdataclass (*CryLasAttenuatorConfig attribute*), 99
is_configdataclass (*CryLasLaserConfig attribute*), 104
is_configdataclass (*EmptyConfig attribute*), 97, 191
is_configdataclass (*FuGConfig attribute*), 113
is_configdataclass (*HeinzingerConfig attribute*), 123
is_configdataclass (*ILS2TConfig attribute*), 176
is_configdataclass (*LaudaProRp245eConfig attribute*), 139
is_configdataclass (*LJMCommunicationConfig attribute*), 17
is_configdataclass (*LuminoxConfig attribute*), 180
is_configdataclass (*MBW973Config attribute*), 142
is_configdataclass (*ModbusTcpCommunicationConfig attribute*), 19
is_configdataclass (*NewportSMC100PPConfig attribute*), 153
is_configdataclass (*OpcUaCommunicationConfig attribute*), 21
is_configdataclass (*PfeifferTPGConfig attribute*), 161
is_configdataclass (*SupercubeConfiguration attribute*), 41, 65
is_configdataclass (*T560Config attribute*), 36
is_configdataclass (*TcpCommunicationConfig attribute*), 27
is_configdataclass (*TechnixConfig attribute*), 81
is_configdataclass (*TiePieDeviceConfig attribute*), 83

is_configdataclass (*VisaCommunicationConfig attribute*), 31
is_data_ready_polling_interval_sec (*TiePieDeviceConfig attribute*), 83
is_done() (*MBW973 method*), 141
is_error() (*ExperimentManager method*), 193
is_finished() (*ExperimentManager method*), 193
is_generic_type_hint() (*in module hvl_ccb.utils.typing*), 189
is_in_range() (*ILS2TRegDatatype method*), 178
is_inactive (*CryLasLaser.LaserStatus property*), 101
is_measurement_data_ready() (*TiePieOscilloscope method*), 92
is_open (*LJMCommunication property*), 15
is_open (*OpcUaCommunication property*), 20
is_open (*SerialCommunication property*), 23
is_open (*TelnetCommunication property*), 28
is_polling() (*Poller method*), 182
is_ready (*CryLasLaser.LaserStatus property*), 101
is_running() (*ExperimentManager method*), 193
is_started (*Technix property*), 80
is_triggered() (*TiePieOscilloscope method*), 92
is_valid_scale_range_reversed_str() (*PfeifferTPGConfig.Model method*), 161
keys() (*CryLasLaserSerialCommunicationConfig class method*), 106
keys() (*EmptyConfig class method*), 97, 191
keys() (*FuGConfig class method*), 113
keys() (*FuGSerialCommunicationConfig class method*), 121
keys() (*HeinzingerConfig class method*), 123
keys() (*HeinzingerSerialCommunicationConfig class method*), 127
keys() (*ILS2TConfig class method*), 176
keys() (*ILS2TModbusTcpCommunicationConfig class method*), 177
keys() (*LaudaProRp245eConfig class method*), 139
keys() (*LaudaProRp245eTcpCommunicationConfig class method*), 140
keys() (*LJMCommunicationConfig class method*), 17
keys() (*LuminoxConfig class method*), 180
keys() (*LuminoxSerialCommunicationConfig class method*), 182
keys() (*MBW973Config class method*), 142
keys() (*MBW973SerialCommunicationConfig class method*), 144
keys() (*ModbusTcpCommunicationConfig class method*), 19
keys() (*NewportSMC100PPConfig class method*), 153
keys() (*NewportSMC100PPSerialCommunicationConfig class method*), 156
keys() (*OpcUaCommunicationConfig class method*), 21
keys() (*PfeifferTPGConfig class method*), 161
keys() (*PfeifferTPGSerialCommunicationConfig class method*), 163
keys() (*PSI9000Config class method*), 109
keys() (*PSI9000VisaCommunicationConfig class method*), 111
keys() (*RTO1024Config class method*), 170
keys() (*RTO1024VisaCommunicationConfig class method*), 171
keys() (*SerialCommunicationConfig class method*), 24
keys() (*SupercubeAOpcUaConfiguration class method*), 58, 76
keys() (*SupercubeBOpcUaConfiguration class method*), 61
keys() (*SupercubeConfiguration class method*), 41, 65
keys() (*SupercubeOpcUaCommunicationConfig class method*), 42, 66
keys() (*T560CommunicationConfig class method*), 34
keys() (*T560Config class method*), 36
keys() (*TcpCommunicationConfig class method*), 27
keys() (*TechnixConfig class method*), 81
keys() (*TechnixSerialCommunicationConfig class method*), 78
keys() (*TechnixTelnetCommunicationConfig class method*), 79
keys() (*TelnetCommunicationConfig class method*), 28

J

J (*LabJack.ThermocoupleType attribute*), 130
jerk_time (*NewportSMC100PPConfig attribute*), 153
JOG (*ILS2T.Mode attribute*), 172
jog_run() (*ILS2T method*), 174
jog_stop() (*ILS2T method*), 174
JOGGING (*NewportStates attribute*), 157
JOGGING_FROM_DISABLE
portSMC100PP.StateMessages
146
JOGGING_FROM_READY
portSMC100PP.StateMessages
146
JOGN_FAST (*ILS2TRegAddr attribute*), 177
JOGN_SLOW (*ILS2TRegAddr attribute*), 177
JR (*NewportConfigCommands attribute*), 145

K

K (*LabJack.TemperatureUnit attribute*), 130
K (*LabJack.ThermocoupleType attribute*), 130
K (*Temperature attribute*), 187
KELVIN (*Temperature attribute*), 187
keys() (*AsyncCommunicationProtocolConfig class method*), 13
keys() (*CryLasAttenuatorConfig class method*), 99
keys() (*CryLasAttenuatorSerialCommunicationConfig class method*), 100
keys() (*CryLasLaserConfig class method*), 104

keys() (*CryLasLaserSerialCommunicationConfig class method*), 106
keys() (*EmptyConfig class method*), 97, 191
keys() (*FuGConfig class method*), 113
keys() (*FuGSerialCommunicationConfig class method*), 121
keys() (*HeinzingerConfig class method*), 123
keys() (*HeinzingerSerialCommunicationConfig class method*), 127
keys() (*ILS2TConfig class method*), 176
keys() (*ILS2TModbusTcpCommunicationConfig class method*), 177
keys() (*LaudaProRp245eConfig class method*), 139
keys() (*LaudaProRp245eTcpCommunicationConfig class method*), 140
keys() (*LJMCommunicationConfig class method*), 17
keys() (*LuminoxConfig class method*), 180
keys() (*LuminoxSerialCommunicationConfig class method*), 182
keys() (*MBW973Config class method*), 142
keys() (*MBW973SerialCommunicationConfig class method*), 144
keys() (*ModbusTcpCommunicationConfig class method*), 19
keys() (*NewportSMC100PPConfig class method*), 153
keys() (*NewportSMC100PPSerialCommunicationConfig class method*), 156
keys() (*OpcUaCommunicationConfig class method*), 21
keys() (*PfeifferTPGConfig class method*), 161
keys() (*PfeifferTPGSerialCommunicationConfig class method*), 163
keys() (*PSI9000Config class method*), 109
keys() (*PSI9000VisaCommunicationConfig class method*), 111
keys() (*RTO1024Config class method*), 170
keys() (*RTO1024VisaCommunicationConfig class method*), 171
keys() (*SerialCommunicationConfig class method*), 24
keys() (*SupercubeAOpcUaConfiguration class method*), 58, 76
keys() (*SupercubeBOpcUaConfiguration class method*), 61
keys() (*SupercubeConfiguration class method*), 41, 65
keys() (*SupercubeOpcUaCommunicationConfig class method*), 42, 66
keys() (*T560CommunicationConfig class method*), 34
keys() (*T560Config class method*), 36
keys() (*TcpCommunicationConfig class method*), 27
keys() (*TechnixConfig class method*), 81
keys() (*TechnixSerialCommunicationConfig class method*), 78
keys() (*TechnixTelnetCommunicationConfig class method*), 79
keys() (*TelnetCommunicationConfig class method*), 28

keys() (*TiePieDeviceConfig class method*), 83
 keys() (*VisaCommunicationConfig class method*), 31
 keys() (*VisaDeviceConfig class method*), 185
 kV (*HeinzingerPNC.UnitVoltage attribute*), 126

L

LabJack (*class in hvl_ccb.dev.labjack*), 128
 LabJack.AInRange (*class in hvl_ccb.dev.labjack*), 129
 LabJack.BitLimit (*class in hvl_ccb.dev.labjack*), 129
 LabJack.CalMicroAmpere (*class in hvl_ccb.dev.labjack*), 129
 LabJack.CjcType (*class in hvl_ccb.dev.labjack*), 129
 LabJack.ClockFrequency (*class in hvl_ccb.dev.labjack*), 129
 LabJack.DeviceType (*class in hvl_ccb.dev.labjack*), 130
 LabJack.DIOStatus (*class in hvl_ccb.dev.labjack*), 129
 LabJack.TemperatureUnit (*class in hvl_ccb.dev.labjack*), 130
 LabJack.ThermocoupleType (*class in hvl_ccb.dev.labjack*), 130
 LabJackError, 134
 LabJackIdentifierDIOError, 134
 laser_off() (*CryLasLaser method*), 101
 laser_on() (*CryLasLaser method*), 102
 LaudaProRp245e (*class in hvl_ccb.dev.lauda*), 134
 LaudaProRp245eCommand (*class in hvl_ccb.dev.lauda*), 136
 LaudaProRp245eCommandError, 138
 LaudaProRp245eConfig (*class in hvl_ccb.dev.lauda*), 138
 LaudaProRp245eConfig.ExtControlModeEnum (*class in hvl_ccb.dev.lauda*), 138
 LaudaProRp245eConfig.OperationModeEnum (*class in hvl_ccb.dev.lauda*), 138
 LaudaProRp245eTcpCommunication (*class in hvl_ccb.dev.lauda*), 139
 LaudaProRp245eTcpCommunicationConfig (*class in hvl_ccb.dev.lauda*), 140
 LEM4000S (*class in hvl_ccb.utils.conversion_sensor*), 185
 LF (*FuGTerminators attribute*), 122
 LFCR (*FuGTerminators attribute*), 122
 line_1 (*MessageBoard attribute*), 54
 line_10 (*MessageBoard attribute*), 54
 line_11 (*MessageBoard attribute*), 54
 line_12 (*MessageBoard attribute*), 54
 line_13 (*MessageBoard attribute*), 54
 line_14 (*MessageBoard attribute*), 54
 line_15 (*MessageBoard attribute*), 54
 line_2 (*MessageBoard attribute*), 54
 line_3 (*MessageBoard attribute*), 54
 line_4 (*MessageBoard attribute*), 54
 line_5 (*MessageBoard attribute*), 54
 line_6 (*MessageBoard attribute*), 54

line_7 (*MessageBoard attribute*), 54
 line_8 (*MessageBoard attribute*), 55
 line_9 (*MessageBoard attribute*), 55
 list_devices() (*TiePieOscilloscope static method*), 92
 list_directory() (*RTO1024 method*), 166
 live (*OpcControl attribute*), 55

LJMCommunication (*class in hvl_ccb.comm.labjack_ljm*), 15
 LJMCommunicationConfig (*class in hvl_ccb.comm.labjack_ljm*), 16
 LJMCommunicationConfig.ConnectionType (*class in hvl_ccb.comm.labjack_ljm*), 16
 LJMCommunicationConfig.DeviceType (*class in hvl_ccb.comm.labjack_ljm*), 16
 LJMCommunicationError, 17
 lm34 (*LabJack.CjcType attribute*), 129
 LMT70A (*class in hvl_ccb.utils.conversion_sensor*), 186
 load_configuration() (*RTO1024 method*), 167
 load_device_configuration() (*T560 method*), 36
 local_display() (*RTO1024 method*), 167
 locked (*DoorStatus attribute*), 49, 68
 LOW (*LabJack.DIOStatus attribute*), 130
 LOWER_TEMP (*LaudaProRp245eCommand attribute*), 137
 lower_temp (*LaudaProRp245eConfig attribute*), 139
 Luminox (*class in hvl_ccb.dev.sst_luminox*), 178
 LuminoxConfig (*class in hvl_ccb.dev.sst_luminox*), 179
 LuminoxException, 180
 LuminoxMeasurementType (*class in hvl_ccb.dev.sst_luminox*), 180
 LuminoxMeasurementTypeDict (*in module hvl_ccb.dev.sst_luminox*), 181
 LuminoxMeasurementTypeError, 181
 LuminoxMeasurementTypeValue (*in module hvl_ccb.dev.sst_luminox*), 181
 LuminoxOutputMode (*class in hvl_ccb.dev.sst_luminox*), 181
 LuminoxOutputModeError, 181
 LuminoxSerialCommunication (*class in hvl_ccb.dev.sst_luminox*), 181
 LuminoxSerialCommunicationConfig (*class in hvl_ccb.dev.sst_luminox*), 181
 LUT (*LMT70A attribute*), 186

M

mA (*HeinzingerPNC.UnitCurrent attribute*), 125
 manual (*EarthingStickOperatingStatus attribute*), 51
 manual() (*EarthingStick class method*), 50, 68
 manual_1 (*EarthingStick attribute*), 50, 68
 manual_2 (*EarthingStick attribute*), 50, 68
 manual_3 (*EarthingStick attribute*), 50, 68
 manual_4 (*EarthingStick attribute*), 50, 68
 manual_5 (*EarthingStick attribute*), 50, 68
 manual_6 (*EarthingStick attribute*), 50, 68
 manuals() (*EarthingStick class method*), 50

<code>MARK (SerialCommunicationParity attribute)</code> , 25		
<code>max_current (FuG property)</code> , 112		
<code>max_current (HeinzingerPNC property)</code> , 126		
<code>max_current (Technix property)</code> , 80		
<code>max_current (TechnixConfig attribute)</code> , 82		
<code>max_current.hardware (FuG property)</code> , 112		
<code>max_current.hardware (HeinzingerPNC property)</code> , 126		
<code>max_pr_number (LaudaProRp245eConfig attribute)</code> , 139		
<code>max_pump_level (LaudaProRp245eConfig attribute)</code> , 139		
<code>max_timeout_retry_nr (OpcUaCommunicationConfig attribute)</code> , 21		
<code>max_voltage (FuG property)</code> , 112		
<code>max_voltage (HeinzingerPNC property)</code> , 126		
<code>max_voltage (Technix property)</code> , 80		
<code>max_voltage (TechnixConfig attribute)</code> , 82		
<code>max_voltage.hardware (FuG property)</code> , 112		
<code>max_voltage.hardware (HeinzingerPNC property)</code> , 126		
<code>MAXIMUM (LabJack.ClockFrequency attribute)</code> , 129		
<code>mbar (PfeifferTPG.PressureUnits attribute)</code> , 158		
<code>MBW973 (class in hvl_ccb.dev.mbw973)</code> , 141		
<code>MBW973Config (class in hvl_ccb.dev.mbw973)</code> , 142		
<code>MBW973ControlRunningException</code> , 143		
<code>MBW973Error</code> , 143		
<code>MBW973PumpRunningException</code> , 143		
<code>MBW973SerialCommunication (class in hvl_ccb.dev.mbw973)</code> , 143		
<code>MBW973SerialCommunicationConfig (class in hvl_ccb.dev.mbw973)</code> , 143		
<code>measure() (PfeifferTPG method)</code> , 160		
<code>measure_all() (PfeifferTPG method)</code> , 160		
<code>measure_current() (HeinzingerDI method)</code> , 124		
<code>measure_voltage() (HeinzingerDI method)</code> , 124		
<code>measure_voltage_current() (PSI9000 method)</code> , 108		
<code>MeasurementsDividerRatio (class in hvl_ccb.dev.supercube.constants)</code> , 54		
<code>MeasurementsDividerRatio (class in hvl_ccb.dev.supercube2015.constants)</code> , 71		
<code>MeasurementsScaledInput (class in hvl_ccb.dev.supercube.constants)</code> , 54		
<code>MeasurementsScaledInput (class in hvl_ccb.dev.supercube2015.constants)</code> , 71		
<code>message (Errors attribute)</code> , 52		
<code>MessageBoard (class in hvl_ccb.dev.supercube.constants)</code> , 54		
<code>micro_step_per_full_step_factor (NewportSMC100PPConfig attribute)</code> , 153		
<code>Micron (PfeifferTPG.PressureUnits attribute)</code> , 158		
<code>MILLIMETER_MERCURY (Pressure attribute)</code> , 187		
<code>MINIMUM (LabJack.ClockFrequency attribute)</code> , 129		
<code>MMHG (Pressure attribute)</code> , 187		
<code>ModbusTcpCommunication (class in hvl_ccb.comm.modbus_tcp)</code> , 18		
<code>ModbusTcpCommunicationConfig (class in hvl_ccb.comm.modbus_tcp)</code> , 19		
<code>ModbusTcpConnectionFailedException</code> , 19		
<code>model (PfeifferTPGConfig attribute)</code> , 161		
<code>module</code>		
<code>hvl_ccb</code> , 194		
<code>hvl_ccb.comm</code> , 32		
<code>hvl_ccb.comm.base</code> , 11		
<code>hvl_ccb.comm.labjack_ljm</code> , 15		
<code>hvl_ccb.comm.modbus_tcp</code> , 18		
<code>hvl_ccb.comm.opc</code> , 20		
<code>hvl_ccb.comm.serial</code> , 22		
<code>hvl_ccb.comm.tcp</code> , 26		
<code>hvl_ccb.comm.telnet</code> , 27		
<code>hvl_ccb.comm.visa</code> , 29		
<code>hvl_ccb.configuration</code> , 190		
<code>hvl_ccb.dev</code> , 185		
<code>hvl_ccb.dev.base</code> , 95		
<code>hvl_ccb.dev.crylas</code> , 97		
<code>hvl_ccb.dev.ea_psi9000</code> , 107		
<code>hvl_ccb.dev.fug</code> , 111		
<code>hvl_ccb.dev.heinzinger</code> , 122		
<code>hvl_ccb.dev.highland_t560</code> , 37		
<code>hvl_ccb.dev.highland_t560.base</code> , 32		
<code>hvl_ccb.dev.highland_t560.channel</code> , 34		
<code>hvl_ccb.dev.highland_t560.device</code> , 35		
<code>hvl_ccb.dev.labjack</code> , 128		
<code>hvl_ccb.dev.lauda</code> , 134		
<code>hvl_ccb.dev.mbw973</code> , 141		
<code>hvl_ccb.dev.newport</code> , 144		
<code>hvl_ccb.dev.pfeiffer_tpg</code> , 158		
<code>hvl_ccb.dev.rs_rto1024</code> , 164		
<code>hvl_ccb.dev.se_ilst2t</code> , 172		
<code>hvl_ccb.dev.sst_luminox</code> , 178		
<code>hvl_ccb.dev.supercube</code> , 61		
<code>hvl_ccb.dev.supercube.base</code> , 37		
<code>hvl_ccb.dev.supercube.constants</code> , 43		
<code>hvl_ccb.dev.supercube.typ_a</code> , 57		
<code>hvl_ccb.dev.supercube.typ_b</code> , 60		
<code>hvl_ccb.dev.supercube2015</code> , 76		
<code>hvl_ccb.dev.supercube2015.base</code> , 61		
<code>hvl_ccb.dev.supercube2015.constants</code> , 67		
<code>hvl_ccb.dev.supercube2015.typ_a</code> , 74		
<code>hvl_ccb.dev.technix</code> , 82		
<code>hvl_ccb.dev.technix.base</code> , 77		
<code>hvl_ccb.dev.technix.device</code> , 80		
<code>hvl_ccb.dev.tiepie</code> , 94		
<code>hvl_ccb.dev.tiepie.base</code> , 83		
<code>hvl_ccb.dev.tiepie.channel</code> , 85		
<code>hvl_ccb.dev.tiepie.device</code> , 87		
<code>hvl_ccb.dev.tiepie.generator</code> , 88		
<code>hvl_ccb.dev.tiepie.i2c</code> , 90		

<code>hvl_ccb.dev.tiepie.oscilloscope</code> , 91	<code>NewportConfigCommands</code>	(class	<i>in</i>
<code>hvl_ccb.dev.tiepie.utils</code> , 93	<i>hvl_ccb.dev.newport)</i> , 144		
<code>hvl_ccb.dev.utils</code> , 182	<code>NewportControllerError</code> , 145		
<code>hvl_ccb.dev.visa</code> , 183	<code>NewportError</code> , 145		
<code>hvl_ccb.exception</code> , 192	<code>NewportMotorError</code> , 145		
<code>hvl_ccb.experiment_manager</code> , 192	<code>NewportMotorPowerSupplyWasCutError</code> , 145		
<code>hvl_ccb.utils</code> , 190	<code>NewportSerialCommunicationError</code> , 157		
<code>hvl_ccb.utils.conversion_sensor</code> , 185	<code>NewportSMC100PP</code> (class <i>in hvl_ccb.dev.newport</i>), 145		
<code>hvl_ccb.utils.conversion_unit</code> , 186	<code>NewportSMC100PP.MotorErrors</code>	(class	<i>in</i>
<code>hvl_ccb.utils.enum</code> , 188	<i>hvl_ccb.dev.newport)</i> , 145		
<code>hvl_ccb.utils.typing</code> , 189	<code>NewportSMC100PP.StateMessages</code>	(class	<i>in</i>
<code>hvl_ccb.utils.validation</code> , 189	<i>hvl_ccb.dev.newport), 146</i>		
<code>MONITOR_I</code> (<i>FuGProbusVRegisterGroups</i> attribute), 119	<code>NewportSMC100PPConfig</code>	(class	<i>in</i>
<code>MONITOR_V</code> (<i>FuGProbusVRegisterGroups</i> attribute), 119	<i>hvl_ccb.dev.newport), 151</i>		
<code>most_recent_error</code> (<i>FuGProbusVConfigRegisters</i> property), 117	<code>NewportSMC100PPConfig.EspStageConfig</code> (class <i>in</i>		
<code>motion_distance_per_full_step</code> (<i>NewportSMC100PPConfig</i> attribute), 153	<i>hvl_ccb.dev.newport), 152</i>		
<code>motor_config</code> (<i>NewportSMC100PPConfig</i> property), 153	<code>NewportSMC100PPSerialCommunication</code> (class <i>in</i>		
<code>move_to_absolute_position()</code> (<i>NewportSMC100PP</i> method), 149	<i>hvl_ccb.dev.newport), 153</i>		
<code>move_to_relative_position()</code> (<i>NewportSMC100PP</i> method), 149	<code>NewportSMC100PPSerialCommunication.ControllerErrors</code> (class <i>in</i>		
<code>move_wait_sec</code> (<i>NewportSMC100PPConfig</i> attribute), 153	<i>hvl_ccb.dev.newport), 154</i>		
<code>MOVING</code> (<i>NewportSMC100PP.StateMessages</i> attribute), 146	<code>NewportSMC100PPSerialCommunicationConfig</code> (class <i>in</i>		
<code>MOVING</code> (<i>NewportStates</i> attribute), 157	<i>hvl_ccb.dev.newport), 156</i>		
<code>MS_NOMINAL_CURRENT</code> (<i>PSI9000</i> attribute), 107	<code>NewportStates</code> (class <i>in hvl_ccb.dev.newport</i>), 157		
<code>MS_NOMINAL_VOLTAGE</code> (<i>PSI9000</i> attribute), 107	<code>NewportUncertainPositionError</code> , 157		
<code>MULTI_COMMANDS_MAX</code> (<i>VisaCommunication</i> attribute), 29	<code>NO</code> (<i>FuDigitalVal</i> attribute), 114		
<code>MULTI_COMMANDS_SEPARATOR</code> (<i>VisaCommunication</i> attribute), 29	<code>NO_ERROR</code> (<i>NewportSMC100PPSerialCommunication.ControllerErrors</i> attribute), 154		
N	<code>NO_REF</code> (<i>NewportStates</i> attribute), 157		
<code>n_channels</code> (<i>TiePieOscilloscope</i> property), 92	<code>NO_REF_ESP_STAGE_ERROR</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
<code>n_max_try_get_device</code> (<i>TiePieDeviceConfig</i> attribute), 83	<i>146</i>		
<code>NameEnum</code> (class <i>in hvl_ccb.utils.enum</i>), 188	<code>NO_REF_FROM_CONFIG</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
<code>NAMES</code> (<i>SerialCommunicationParity</i> attribute), 25	<i>146</i>		
<code>names()</code> (<i>RTO1024.TriggerModes</i> class method), 165	<code>NO_REF_FROM_DISABLED</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
<code>namespace_index</code> (<i>SupercubeConfiguration</i> attribute), 42, 65	<i>146</i>		
<code>NATIVEONLY</code> (<i>TiePieOscilloscopeAutoResolutionModes</i> attribute), 92	<code>NO_REF_FROM_HOMING</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
<code>NED_END_OF_TURN</code> (<i>NewportSMC100PP.MotorErrors</i> attribute), 145	<i>146</i>		
<code>NEG</code> (<i>ILS2T.Refl6Jog</i> attribute), 172	<code>NO_REF_FROM_JOGGING</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
<code>NEG_FAST</code> (<i>ILS2T.Refl6Jog</i> attribute), 172	<i>146</i>		
<code>NEGATIVE</code> (<i>FuGPolarities</i> attribute), 116	<code>NO_REF_FROM_MOVING</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
<code>negative_software_limit</code> (<i>NewportSMC100PPConfig</i> attribute), 153	<i>146</i>		
	<code>NO_REF_FROM_READY</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
	<i>146</i>		
	<code>NO_REF_FROM_RESET</code>	(<i>NewportSMC100PP.StateMessages</i> attribute),	
	<i>146</i>		

No_sensor (*PfeifferTPG.SensorStatus attribute*), 158
 NOISE (*TiePieGeneratorSignalType attribute*), 89
 NONE (*ILS2T.Ref16Jog attribute*), 172
 NONE (*LabJack.ThermocoupleType attribute*), 130
 None (*PfeifferTPG.SensorTypes attribute*), 159
 NONE (*SerialCommunicationParity attribute*), 25
 NoPower (*PowerSetup attribute*), 56
 NORMAL (*RTO1024.TriggerModes attribute*), 165
 noSen (*PfeifferTPG.SensorTypes attribute*), 159
 noSENSOR (*PfeifferTPG.SensorTypes attribute*), 159
 not_defined (*AlarmText attribute*), 44, 67
 nr_trials_activate (*LuminoxConfig attribute*), 180
 NullCommunicationProtocol (class in *hvl_ccb.comm.base*), 14
 number (*EarthingStick property*), 50
 Number (in module *hvl_ccb.utils.typing*), 189
 number_of_decimals (*HeinzingerConfig attribute*), 123
 number_of_sensors (*PfeifferTPG property*), 160

O

ODD (*SerialCommunicationParity attribute*), 25
 OFF (*AutoInstallMode attribute*), 33
 OFF (*FuGDigitalVal attribute*), 114
 OFF (*GateMode attribute*), 33
 OFF (*HeinzingerDI.OutputStatus attribute*), 124
 OFF (*TriggerMode attribute*), 34
 offset (*TiePieGeneratorConfig property*), 89
 OH (*NewportConfigCommands attribute*), 145
 Ok (*PfeifferTPG.SensorStatus attribute*), 159
 on (*FuG property*), 112
 ON (*FuGDigitalVal attribute*), 114
 on (*FuGProbusVDIRegisters property*), 118
 ON (*HeinzingerDI.OutputStatus attribute*), 124
 ON (*ILS2T.State attribute*), 173
 ONE (*HeinzingerConfig.RecordingsEnum attribute*), 123
 ONE (*LabJack.AInRange attribute*), 129
 ONE (*SerialCommunicationStopbits attribute*), 25
 ONE_HUNDREDTH (*LabJack.AInRange attribute*), 129
 ONE_POINT_FIVE (*SerialCommunicationStopbits attribute*), 25
 ONE_TENTH (*LabJack.AInRange attribute*), 129
 ONLYUPWARDSOFFTOZERO (*FuGRampModes attribute*), 120
 OpcControl (class in *hvl_ccb.dev.supercube.constants*), 55
 OpcUaCommunication (class in *hvl_ccb.comm.opc*), 20
 OpcUaCommunicationConfig (class in *hvl_ccb.comm.opc*), 21
 OpcUaCommunicationIOError, 22
 OpcUaCommunicationTimeoutError, 22
 OpcUaSubHandler (class in *hvl_ccb.comm.opc*), 22
 open (*DoorStatus attribute*), 49, 68
 open (*EarthingStickOperation attribute*), 51
 open (*EarthingStickStatus attribute*), 52, 70

open() (*CommunicationProtocol method*), 14
 open() (*LaudaProRp245eTcpCommunication method*), 139
 open() (*LJMCommunication method*), 15
 open() (*ModbusTcpCommunication method*), 18
 open() (*NullCommunicationProtocol method*), 14
 open() (*OpcUaCommunication method*), 20
 open() (*SerialCommunication method*), 23
 open() (*Tcp method*), 26
 open() (*TelnetCommunication method*), 28
 open() (*VisaCommunication method*), 30
 open_interlock (*Technix property*), 81
 open_shutter() (*CryLasLaser method*), 102
 open_timeout (*VisaCommunicationConfig attribute*), 31
 OPENED (*CryLasLaser.AnswersShutter attribute*), 100
 OPENED (*CryLasLaserShutterStatus attribute*), 106
 operate() (*Supercube2015Base method*), 63
 operate() (*SupercubeBase method*), 39
 operate_earthing_stick() (*SupercubeBase method*), 39
 operating_status() (*EarthingStick class method*), 50
 operating_status_1 (*EarthingStick attribute*), 50
 operating_status_2 (*EarthingStick attribute*), 50
 operating_status_3 (*EarthingStick attribute*), 50
 operating_status_4 (*EarthingStick attribute*), 50
 operating_status_5 (*EarthingStick attribute*), 51
 operating_status_6 (*EarthingStick attribute*), 51
 operating_statuses() (*EarthingStick class method*), 51
 OPERATION_MODE (*LaudaProRp245eCommand attribute*), 137
 operation_mode (*LaudaProRp245eConfig attribute*), 139
 optional_defaults() (*AsyncCommunicationProtocolConfig class method*), 13
 optional_defaults() (*CryLasAttenuatorConfig class method*), 99
 optional_defaults() (*CryLasAttenuatorSerialCommunicationConfig class method*), 100
 optional_defaults() (*CryLasLaserConfig class method*), 104
 optional_defaults() (*CryLasLaserSerialCommunicationConfig class method*), 106
 optional_defaults() (*EmptyConfig class method*), 97, 191
 optional_defaults() (*FuGConfig class method*), 113
 optional_defaults() (*FuGSerialCommunicationConfig class method*), 121
 optional_defaults() (*HeinzingerConfig class method*), 123
 optional_defaults() (*HeinzingerSerialCommunicationConfig class method*), 127
 optional_defaults() (*ILS2TConfig class method*), 176

optional_defaults()
 (ILS2TModbusTcpCommunicationConfig
 class method), 177

optional_defaults() (LaudaProRp245eConfig class
 method), 139

optional_defaults() (LaudaProRp245eTcpCommunicationConfig class
 method), 140

optional_defaults() (LJMCommunicationConfig
 class method), 17

optional_defaults() (LuminoxConfig class method),
 180

optional_defaults() (LuminoxSerialCommunicationConfig class method), 182

optional_defaults() (MBW973Config class method),
 142

optional_defaults()
 (MBW973SerialCommunicationConfig class
 method), 144

optional_defaults() (ModbusTcpCommunicationConfig class method), 19

optional_defaults() (NewportSMC100PPConfig class method), 153

optional_defaults() (NewportSMC100PPSerialCommunicationConfig
 class method), 157

optional_defaults() (OpcUaCommunicationConfig
 class method), 21

optional_defaults() (PfeifferTPGConfig class
 method), 161

optional_defaults() (PfeifferTPGSerialCommunicationConfig class method), 163

optional_defaults() (PSI9000Config class method),
 109

optional_defaults()
 (PSI9000VisaCommunicationConfig class
 method), 111

optional_defaults() (RTO1024Config class method),
 170

optional_defaults()
 (RTO1024VisaCommunicationConfig class
 method), 171

optional_defaults() (SerialCommunicationConfig
 class method), 24

optional_defaults() (SupercubeAOpcUaConfiguration class method), 58, 76

optional_defaults() (SupercubeBOpcUaConfiguration class method), 61

optional_defaults() (SupercubeConfiguration class
 method), 42, 65

optional_defaults() (SupercubeOpcUaCommunicationConfig class method), 42, 66

optional_defaults() (T560CommunicationConfig class method), 34

optional_defaults() (T560Config class method), 36

optional_defaults() (TcpCommunicationConfig
 class method), 27

optional_defaults() (TechnixConfig class method),
 82

optional_defaults() (TechnixSerialCommunicationConfig class method), 78

optional_defaults() (TechnixTelnetCommunicationConfig class method), 79

optional_defaults() (TelnetCommunicationConfig
 class method), 28

optional_defaults() (TiePieDeviceConfig class
 method), 84

optional_defaults() (VisaCommunicationConfig
 class method), 31

optional_defaults() (VisaDeviceConfig class
 method), 185

OSCILLOSCOPE (TiePieDeviceType attribute), 84

OT (NewportConfigCommands attribute), 145

out (FuGProbusVDORegisters property), 118

out_1_1 (GeneralSupport attribute), 53, 71

out_1_2 (GeneralSupport attribute), 53, 71

out_2_1 (GeneralSupport attribute), 53, 71

out_2_2 (GeneralSupport attribute), 53, 71

out_3_1 (GeneralSupport attribute), 53, 71

out_3_2 (GeneralSupport attribute), 53, 71

out_4_1 (GeneralSupport attribute), 53, 71

out_4_2 (GeneralSupport attribute), 53, 71

out_5_1 (GeneralSupport attribute), 53, 71

out_5_2 (GeneralSupport attribute), 53, 71

out_6_1 (GeneralSupport attribute), 53, 71

out_6_2 (GeneralSupport attribute), 53, 71

OUTPUT (FuGProbusIVCommands attribute), 117

OUTPUT (GateMode attribute), 33

output (Technix property), 81

output() (GeneralSupport class method), 53, 71

output_off() (FuGProbusIV method), 116

output_off() (HeinzingerDI method), 124

output_on() (HeinzingerDI method), 124

OUTPUT_POWER_EXCEEDED (NewportSMC100PP.MotorErrors attribute),
 145

output_status (HeinzingerDI property), 125

OUTPUTONCMD (FuGProbusVRegisterGroups attribute),
 119

OUTPUTX0 (FuGProbusVRegisterGroups attribute), 119

OUTPUTX1 (FuGProbusVRegisterGroups attribute), 119

OUTPUTX2 (FuGProbusVRegisterGroups attribute), 119

OUTPUTXCMD (FuGProbusVRegisterGroups attribute),
 119

outX0 (FuG property), 112

outX1 (FuG property), 112

outX2 (FuG property), 112

outXCMD (FuG property), 112

Overrange (*PfeifferTPG.SensorStatus* attribute), 159

P

PA (*Pressure* attribute), 187

PARAM_MISSING_OR_INVALID (New-
portSMC100PPSerialCommunication.ControllerErrors
attribute), 154

parity (*CryLasAttenuatorSerialCommunicationConfig*
attribute), 100

parity (*CryLasLaserSerialCommunicationConfig*
attribute), 106

parity (*FuGSerialCommunicationConfig* attribute), 121

parity (*HeinzingerSerialCommunicationConfig* at-
tribute), 127

parity (*LuminoxSerialCommunicationConfig* attribute),
182

parity (*MBW973SerialCommunicationConfig* at-
tribute), 144

parity (*NewportSMC100PPSerialCommunicationConfig*
attribute), 157

parity (*PfeifferTPGSerialCommunicationConfig* at-
tribute), 163

Parity (*SerialCommunicationConfig* attribute), 24

parity (*SerialCommunicationConfig* attribute), 24

parse_read_measurement_value() (*LuminoxMea-
surementType* method), 180

partial_pressure_o2 (*LuminoxMeasurementType* at-
tribute), 180

Pascal (*PfeifferTPG.PressureUnits* attribute), 158

PASCAL (*Pressure* attribute), 187

pause() (*LaudaProRp245e* method), 135

pause_ramp() (*LaudaProRp245e* method), 135

PBR (*PfeifferTPG.SensorTypes* attribute), 159

PEAK_CURRENT_LIMIT (New-
portSMC100PP.MotorErrors
attribute), 146

peak_output_current_limit (New-
portSMC100PPConfig attribute), 153

percent_o2 (*LuminoxMeasurementType* attribute), 180

period (*T560* property), 36

PfeifferTPG (class in *hvl_ccb.dev.pfeiffer_tpg*), 158

PfeifferTPG.PressureUnits (class
hvl_ccb.dev.pfeiffer_tpg), 158

PfeifferTPG.SensorStatus (class
hvl_ccb.dev.pfeiffer_tpg), 158

PfeifferTPG.SensorTypes (class
hvl_ccb.dev.pfeiffer_tpg), 159

PfeifferTPGConfig (class in *hvl_ccb.dev.pfeiffer_tpg*),
161

PfeifferTPGConfig.Model (class
hvl_ccb.dev.pfeiffer_tpg), 161

PfeifferTPGError, 162

PfeifferTPGSerialCommunication (class
hvl_ccb.dev.pfeiffer_tpg), 162

PfeifferTPGSerialCommunicationConfig (class in
hvl_ccb.dev.pfeiffer_tpg), 162

PKR (*PfeifferTPG.SensorTypes* attribute), 159

Polarity (class in *hvl_ccb.dev.highland_t560.base*), 33

POLARITY (*FuGProbusIVCommands* attribute), 117

Poller (class in *hvl_ccb.dev.utils*), 182

polling (*LuminoxOutputMode* attribute), 181

polling_delay_sec (*SupercubeConfiguration* at-
tribute), 42

polling_interval (*MBW973Config* attribute), 143

polling_interval_sec (*SupercubeConfiguration* at-
tribute), 42

polling_interval_sec (*TechnixConfig* attribute), 82

polling_period (*CryLasLaserConfig* attribute), 104

polling_timeout (*CryLasLaserConfig* attribute), 104

port (*ModbusTcpCommunicationConfig* attribute), 19

port (*OpcUaCommunicationConfig* attribute), 21

port (*SerialCommunicationConfig* attribute), 24

port (*SupercubeOpcUaCommunicationConfig* attribute),
66

port (*T560CommunicationConfig* attribute), 34

port (*TcpCommunicationConfig* attribute), 27

port (*TechnixTelnetCommunicationConfig* attribute), 79

port (*TelnetCommunicationConfig* attribute), 29

port (*VisaCommunicationConfig* attribute), 31

port_range() (*GeneralSupport* class method), 53

POS (*ILS2T.Ref16Jog* attribute), 173

POS_END_OF_TURN (*NewportSMC100PP.MotorErrors*
attribute), 146

POS_FAST (*ILS2T.Ref16Jog* attribute), 173

POSITION (*ILS2TRegAddr* attribute), 177

POSITION_OUT_OF_LIMIT (New-
portSMC100PPSerialCommunication.ControllerErrors
attribute), 154

POSITIVE (*FuGPolarities* attribute), 116

positive_software_limit (New-
portSMC100PPConfig attribute), 153

post_force_value() (*NewportSMC100PPConfig*
method), 153

post_stop_pause_sec (*TechnixConfig* attribute), 82

POUNDS_PER_SQUARE_INCH (*Pressure* attribute), 187

Power (class in *hvl_ccb.dev.supercube.constants*), 55

Power (class in *hvl_ccb.dev.supercube2015.constants*),
72

power_limit (*PSI9000Config* attribute), 110

PowerSetup (class in *hvl_ccb.dev.supercube.constants*),
55

PowerSetup (class in *hvl_ccb.dev.supercube2015.constants*),
72

pre_sample_ratio (*TiePieOscilloscopeConfig* prop-
erty), 93

prepare_ultra_segmentation() (*RTO1024* method),
167

<p><code>preserve_type()</code> (in module <code>hvl_ccb.utils.conversion_unit</code>), 187</p> <p><code>Pressure</code> (class in <code>hvl_ccb.utils.conversion_unit</code>), 186</p> <p><code>probe_offset</code> (<code>TiePieOscilloscopeChannelConfig</code> property), 86</p> <p><code>PSI</code> (<code>Pressure</code> attribute), 187</p> <p><code>PSI9000</code> (class in <code>hvl_ccb.dev.ea_psi9000</code>), 107</p> <p><code>PSI9000Config</code> (class in <code>hvl_ccb.dev.ea_psi9000</code>), 109</p> <p><code>PSI9000Error</code>, 110</p> <p><code>PSI9000VisaCommunication</code> (class in <code>hvl_ccb.dev.ea_psi9000</code>), 110</p> <p><code>PSI9000VisaCommunicationConfig</code> (class in <code>hvl_ccb.dev.ea_psi9000</code>), 110</p> <p><code>PT100</code> (<code>LabJack.ThermocoupleType</code> attribute), 130</p> <p><code>PT1000</code> (<code>LabJack.ThermocoupleType</code> attribute), 130</p> <p><code>PT500</code> (<code>LabJack.ThermocoupleType</code> attribute), 130</p> <p><code>PTP</code> (<code>ILS2T.Mode</code> attribute), 172</p> <p><code>PublicPropertiesReprMixin</code> (class in <code>hvl_ccb.dev.tiepie.utils</code>), 93</p> <p><code>PULSE</code> (<code>TiePieGeneratorSignalType</code> attribute), 89</p> <p><code>pump_init</code> (<code>LaudaProRp245eConfig</code> attribute), 139</p> <p><code>PUMP_LEVEL</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p>	<p>R</p> <p><code>R</code> (<code>LabJack.ThermocoupleType</code> attribute), 130</p> <p><code>raise_()</code> (<code>FuGErrorcodes</code> method), 115</p> <p><code>RAMP_ACC</code> (<code>ILS2TRegAddr</code> attribute), 177</p> <p><code>RAMP_CONTINUE</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_DECEL</code> (<code>ILS2TRegAddr</code> attribute), 177</p> <p><code>RAMP_DELETE</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_ITERATIONS</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_N_MAX</code> (<code>ILS2TRegAddr</code> attribute), 177</p> <p><code>RAMP_PAUSE</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_SELECT</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_SET</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_START</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_STOP</code> (<code>LaudaProRp245eCommand</code> attribute), 137</p> <p><code>RAMP_TYPE</code> (<code>ILS2TRegAddr</code> attribute), 177</p> <p><code>rampmode</code> (<code>FuGProbusVSetRegisters</code> property), 120</p> <p><code>ramprate</code> (<code>FuGProbusVSetRegisters</code> property), 120</p> <p><code>rampstate</code> (<code>FuGProbusVSetRegisters</code> property), 120</p> <p><code>RAMPUPWARDS</code> (<code>FuGRampModes</code> attribute), 120</p> <p><code>range()</code> (<code>EarthingStick</code> class method), 51</p> <p><code>RATEDCURRENT</code> (<code>FuGReadbackChannels</code> attribute), 120</p> <p><code>RATEDVOLTAGE</code> (<code>FuGReadbackChannels</code> attribute), 120</p> <p><code>read()</code> (<code>AsyncCommunicationProtocol</code> method), 11</p> <p><code>read()</code> (<code>CryLasLaserSerialCommunication</code> method), 105</p> <p><code>read()</code> (<code>LaudaProRp245eTcpCommunication</code> method), 139</p> <p><code>read()</code> (<code>MBW973</code> method), 141</p> <p><code>read()</code> (<code>OpcUaCommunication</code> method), 20</p> <p><code>read()</code> (<code>Supercube2015Base</code> method), 63</p> <p><code>read()</code> (<code>SupercubeBase</code> method), 40</p> <p><code>read()</code> (<code>Tcp</code> method), 26</p> <p><code>read_all()</code> (<code>AsyncCommunicationProtocol</code> method), 11</p> <p><code>read_bytes()</code> (<code>AsyncCommunicationProtocol</code> method), 12</p> <p><code>read_bytes()</code> (<code>SerialCommunication</code> method), 23</p> <p><code>read_bytes()</code> (<code>TelnetCommunication</code> method), 28</p> <p><code>read_float()</code> (<code>MBW973</code> method), 141</p> <p><code>read_holding_registers()</code> (<code>ModbusTcpCommunication</code> method), 18</p> <p><code>read_input_registers()</code> (<code>ModbusTcpCommunication</code> method), 18</p> <p><code>read_int()</code> (<code>MBW973</code> method), 141</p> <p><code>read_measurement()</code> (<code>RTO1024</code> method), 167</p> <p><code>read_measurements()</code> (<code>MBW973</code> method), 142</p> <p><code>read_name()</code> (<code>LJMCommunication</code> method), 15</p> <p><code>read_nonempty()</code> (<code>AsyncCommunicationProtocol</code> method), 12</p> <p><code>read_output_while_polling</code> (<code>TechnixConfig</code> attribute), 82</p>
--	--

read_resistance() (*LabJack method*), 132
read_single_bytes() (*SerialCommunication method*), 23
read_streaming() (*Luminox method*), 179
read_termination (*VisaCommunicationConfig attribute*), 31
read_text() (*AsyncCommunicationProtocol method*), 12
read_text() (*NewportSMC100PPSerialCommunication method*), 155
read_text_nonempty() (*AsyncCommunicationProtocol method*), 12
READ_TEXT_SKIP_PREFIXES (*CryLasLaserSerialCommunication attribute*), 104
read_thermocouple() (*LabJack method*), 132
readback_data (*FuGProbusVConfigRegisters property*), 117
READBACKCHANNEL (*FuGProbusIVCommands attribute*), 117
READY (*CryLasLaser.AnswersStatus attribute*), 101
READY (*ILS2T.State attribute*), 173
READY (*NewportStates attribute*), 157
ready() (*Supercube2015Base method*), 63
ready() (*SupercubeBase method*), 40
READY_ACTIVE (*CryLasLaser.LaserStatus attribute*), 101
READY_FROM_DISABLE (*NewportSMC100PP.StateMessages attribute*), 146
READY_FROM_HOMING (*NewportSMC100PP.StateMessages attribute*), 146
READY_FROM_JOGGING (*NewportSMC100PP.StateMessages attribute*), 146
READY_FROM_MOVING (*NewportSMC100PP.StateMessages attribute*), 146
READY_INACTIVE (*CryLasLaser.LaserStatus attribute*), 101
record_length (*TiePieOscilloscopeConfig property*), 93
RedOperate (*SafetyStatus attribute*), 56, 73
RedReady (*SafetyStatus attribute*), 56, 74
reg_3 (*FuGProbusVDIRegisters property*), 118
RegAddr (*ILS2T attribute*), 173
RegDatatype (*ILS2T attribute*), 173
register_pulse_time (*TechnixConfig attribute*), 82
RELATIVE (*TiePieOscilloscopeTriggerLevelMode attribute*), 87
RELATIVE_POSITION_MOTOR (*ILS2T.ActionsPtp attribute*), 172
RELATIVE_POSITION_TARGET (*ILS2T.ActionsPtp attribute*), 172
remote (*Technix property*), 81
remove_device() (*DeviceSequenceMixin method*), 96
require_block_measurement_support (*TiePieDeviceConfig attribute*), 84
required_keys() (*AsyncCommunicationProtocolConfig class method*), 13
required_keys() (*CryLasAttenuatorConfig class method*), 99
required_keys() (*CryLasAttenuatorSerialCommunicationConfig class method*), 100
required_keys() (*CryLasLaserConfig class method*), 104
required_keys() (*CryLasLaserSerialCommunicationConfig class method*), 106
required_keys() (*EmptyConfig class method*), 97, 191
required_keys() (*FuGConfig class method*), 113
required_keys() (*FuGSerialCommunicationConfig class method*), 121
required_keys() (*HeinzingerConfig class method*), 123
required_keys() (*HeinzingerSerialCommunicationConfig class method*), 127
required_keys() (*ILS2TConfig class method*), 176
required_keys() (*ILS2TModbusTcpCommunicationConfig class method*), 177
required_keys() (*LaudaProRp245eConfig class method*), 139
required_keys() (*LaudaProRp245eTcpCommunicationConfig class method*), 140
required_keys() (*LJMCommunicationConfig class method*), 17
required_keys() (*LuminoxConfig class method*), 180
required_keys() (*LuminoxSerialCommunicationConfig class method*), 182
required_keys() (*MBW973Config class method*), 143
required_keys() (*MBW973SerialCommunicationConfig class method*), 144
required_keys() (*ModbusTcpCommunicationConfig class method*), 19
required_keys() (*NewportSMC100PPConfig class method*), 153
required_keys() (*NewportSMC100PPSerialCommunicationConfig class method*), 157
required_keys() (*OpcUaCommunicationConfig class method*), 22
required_keys() (*PfeifferTPGConfig class method*), 161
required_keys() (*PfeifferTPGSerialCommunicationConfig class method*), 163
required_keys() (*PSI9000Config class method*), 110
required_keys() (*PSI9000VisaCommunicationConfig class method*), 111
required_keys() (*RTO1024Config class method*), 171

<code>required_keys()</code> (<i>RTO1024VisaCommunicationConfig class method</i>), 171	<code>RT01024VisaCommunication</code> (class in <i>hvl_ccb.dev.rs_rto1024</i>), 171
<code>required_keys()</code> (<i>SerialCommunicationConfig class method</i>), 25	<code>RT01024VisaCommunicationConfig</code> (class in <i>hvl_ccb.dev.rs_rto1024</i>), 171
<code>required_keys()</code> (<i>SupercubeAOpcUaConfiguration class method</i>), 58, 76	<code>run()</code> (<i>ExperimentManager method</i>), 193
<code>required_keys()</code> (<i>SupercubeBOpcUaConfiguration class method</i>), 61	<code>run()</code> (<i>LaudaProRp245e method</i>), 135
<code>required_keys()</code> (<i>SupercubeConfiguration class method</i>), 42, 65	<code>run_continuous_acquisition()</code> (<i>RTO1024 method</i>), 167
<code>required_keys()</code> (<i>SupercubeOpcUaCommunicationConfig class method</i>), 43, 66	<code>run_single_acquisition()</code> (<i>RTO1024 method</i>), 167
<code>required_keys()</code> (<i>T560CommunicationConfig class method</i>), 34	<code>RUNNING</code> (<i>ExperimentStatus attribute</i>), 193
<code>required_keys()</code> (<i>T560Config class method</i>), 37	
<code>required_keys()</code> (<i>TcpCommunicationConfig class method</i>), 27	S
<code>required_keys()</code> (<i>TechnixConfig class method</i>), 82	<code>S</code> (<i>LabJack.ThermocoupleType attribute</i>), 131
<code>required_keys()</code> (<i>TechnixSerialCommunicationConfig class method</i>), 78	<code>SA</code> (<i>NewportConfigCommands attribute</i>), 145
<code>required_keys()</code> (<i>TechnixTelnetCommunicationConfig class method</i>), 79	<code>SafeGround</code> (class in <i>hvl_ccb.dev.tiepie.channel</i>), 85
<code>required_keys()</code> (<i>TelnetCommunicationConfig class method</i>), 29	<code>Safety</code> (class in <i>hvl_ccb.dev.supercube.constants</i>), 56
<code>required_keys()</code> (<i>TiePieDeviceConfig class method</i>), 84	<code>Safety</code> (class in <i>hvl_ccb.dev.supercube2015.constants</i>), 73
<code>required_keys()</code> (<i>VisaCommunicationConfig class method</i>), 31	<code>SafetyStatus</code> (class in <i>hvl_ccb.dev.supercube.constants</i>), 56
<code>required_keys()</code> (<i>VisaDeviceConfig class method</i>), 185	<code>SafetyStatus</code> (class in <i>hvl_ccb.dev.supercube2015.constants</i>), 73
<code>reset</code> (<i>BreakdownDetection attribute</i>), 49, 68	<code>sample_frequency</code> (<i>TiePieOscilloscopeConfig property</i>), 93
<code>RESET</code> (<i>FuGProbusIVCommands attribute</i>), 117	<code>save_configuration()</code> (<i>RTO1024 method</i>), 167
<code>reset()</code> (<i>FuGProbusIV method</i>), 116	<code>save_device_configuration()</code> (<i>T560 method</i>), 36
<code>reset()</code> (<i>NewportSMC100PP method</i>), 150	<code>save_waveform_history()</code> (<i>RTO1024 method</i>), 167
<code>reset()</code> (<i>VisaDevice method</i>), 184	<code>SCALE</code> (<i>ILS2TRegAddr attribute</i>), 177
<code>reset_error()</code> (<i>ILS2T method</i>), 174	<code>ScalingFactorValueError</code> , 178
<code>reset_interface()</code> (<i>HeinzingerDI method</i>), 125	<code>screw_scaling</code> (<i>NewportSMC100PPConfig attribute</i>), 153
<code>reset_ramp()</code> (<i>LaudaProRp245e method</i>), 135	<code>send_command()</code> (<i>NewportSMC100PPSerialCommunication method</i>), 155
<code>resolution</code> (<i>TiePieOscilloscopeConfig property</i>), 93	<code>send_command()</code> (<i>PfeifferTPGSerialCommunication method</i>), 162
<code>response_sleep_time</code> (<i>CryLasAttenuatorConfig attribute</i>), 99	<code>send_pulses()</code> (<i>LabJack method</i>), 132
<code>RISING</code> (<i>TiePieOscilloscopeTriggerKind attribute</i>), 87	<code>send_stop()</code> (<i>NewportSMC100PPSerialCommunication method</i>), 156
<code>RISING_OR_FALLING</code> (<i>TiePieOscilloscopeTriggerKind attribute</i>), 87	<code>Sensor</code> (class in <i>hvl_ccb.utils.conversion_sensor</i>), 186
<code>RMS_CURRENT_LIMIT</code> (<i>NewportSMC100PP.MotorErrors attribute</i>), 146	<code>Sensor_error</code> (<i>PfeifferTPG.SensorStatus attribute</i>), 159
<code>rpm_max_init</code> (<i>ILS2TConfig attribute</i>), 176	<code>Sensor_off</code> (<i>PfeifferTPG.SensorStatus attribute</i>), 159
<code>rs485_address</code> (<i>NewportSMC100PPConfig attribute</i>), 153	<code>sensor_status</code> (<i>LuminoxMeasurementType attribute</i>), 180
<code>RT01024</code> (class in <i>hvl_ccb.dev.rs_rto1024</i>), 164	<code>SERIAL</code> (<i>LaudaProRp245eConfig.ExtControlModeEnum attribute</i>), 138
<code>RT01024.TriggerModes</code> (class in <i>hvl_ccb.dev.rs_rto1024</i>), 164	<code>serial_number</code> (<i>LuminoxMeasurementType attribute</i>), 180
<code>RT01024Config</code> (class in <i>hvl_ccb.dev.rs_rto1024</i>), 170	<code>serial_number</code> (<i>TiePieDeviceConfig attribute</i>), 84
<code>RT01024Error</code> , 171	<code>SerialCommunication</code> (class in <i>hvl_ccb.comm.serial</i>), 22
	<code>SerialCommunicationBytesize</code> (class in <i>hvl_ccb.comm.serial</i>), 23

SerialCommunicationConfig	(class	<i>in</i>	set_output() (<i>PSI9000 method</i>), 108
<i>hvl_ccb.comm.serial</i>), 23			set_positive_software_limit() (New- portSMC100PP method), 151
SerialCommunicationIOError	, 25		
SerialCommunicationParity	(class	<i>in</i>	set_pulse_energy() (<i>CryLasLaser method</i>), 102
<i>hvl_ccb.comm.serial</i>), 25			set_pump_level() (<i>LaudaProRp245e method</i>), 135
SerialCommunicationStopbits	(class	<i>in</i>	set_ramp_iterations() (<i>LaudaProRp245e method</i>), 135
<i>hvl_ccb.comm.serial</i>), 25			set_ramp_program() (<i>LaudaProRp245e method</i>), 136
set_acceleration() (<i>NewportSMC100PP method</i>),			set_ramp_segment() (<i>LaudaProRp245e method</i>), 136
150			set_ramp_type() (<i>ILS2T method</i>), 175
set_acquire_length() (<i>RTO1024 method</i>),	167		set_reference_point() (<i>RTO1024 method</i>), 169
set_ain_differential() (<i>LabJack method</i>),	132		set_register() (<i>FuGProbusV method</i>), 117
set_ain_range() (<i>LabJack method</i>),	133		set_remote_control() (<i>Supercube2015Base method</i>), 64
set_ain_resistance() (<i>LabJack method</i>),	133		set_remote_control() (<i>SupercubeBase method</i>), 40
set_ain_resolution() (<i>LabJack method</i>),	133		set_repetition_rate() (<i>CryLasLaser method</i>), 102
set_ain_thermocouple() (<i>LabJack method</i>),	133		set_repetitions() (<i>RTO1024 method</i>), 169
set_analog_output() (<i>LabJack method</i>),	133		set_slope() (<i>Supercube2015WithFU method</i>), 75
set_attenuation() (<i>CryLasAttenuator method</i>),	98		set_slope() (<i>SupercubeWithFU method</i>), 59
set_cee16_socket() (<i>Supercube2015Base method</i>),			set_status_board() (<i>SupercubeBase method</i>), 40
63			set_support_output() (<i>Supercube2015Base method</i>), 64
set_cee16_socket() (<i>SupercubeBase method</i>),	40		set_support_output() (<i>SupercubeBase method</i>), 40
set_channel_offset() (<i>RTO1024 method</i>),	168		set_support_output_impulse() (<i>Supercube2015Base method</i>), 64
set_channel_position() (<i>RTO1024 method</i>),	168		set_support_output_impulse() (<i>SupercubeBase method</i>), 40
set_channel_range() (<i>RTO1024 method</i>),	168		set_system_lock() (<i>PSI9000 method</i>), 108
set_channel_scale() (<i>RTO1024 method</i>),	168		set_t13_socket() (<i>Supercube2015Base method</i>), 64
set_channel_state() (<i>RTO1024 method</i>),	169		set_t13_socket() (<i>SupercubeBase method</i>), 41
set_clock() (<i>LabJack method</i>),	134		set_target_voltage() (<i>Supercube2015WithFU method</i>), 75
set_control_mode() (<i>LaudaProRp245e method</i>),	135		set_target_voltage() (<i>SupercubeWithFU method</i>), 59
set_current() (<i>HeinzingerDI method</i>),	125		set_temp_set_point() (<i>LaudaProRp245e method</i>), 136
set_current() (<i>HeinzingerPNC method</i>),	126		set_transmission() (<i>CryLasAttenuator method</i>), 98
set_digital_output() (<i>LabJack method</i>),	134		set_trigger_level() (<i>RTO1024 method</i>), 169
set_display_unit() (<i>PfeifferTPG method</i>),	160		set_trigger_mode() (<i>RTO1024 method</i>), 170
set_earthing_manual() (<i>Supercube2015Base method</i>),			set_trigger_source() (<i>RTO1024 method</i>), 170
63			set_upper_limits() (<i>PSI9000 method</i>), 108
set_external_temp() (<i>LaudaProRp245e method</i>),	135		set_voltage() (<i>HeinzingerDI method</i>), 125
set_full_scale_mbar() (<i>PfeifferTPG method</i>),	160		set_voltage() (<i>HeinzingerPNC method</i>), 126
set_full_scale_unitless() (<i>PfeifferTPG method</i>),			set_voltage_current() (<i>PSI9000 method</i>), 108
160			SETCURRENT (<i>FuGProbusVRegisterGroups attribute</i>), 119
set_init_attenuation() (<i>CryLasAttenuator method</i>),	98		setup (<i>Power attribute</i>), 55, 72
set_init_shutter_status() (<i>CryLasLaser method</i>),			setvalue (<i>FuGProbusVSetRegisters property</i>), 120
102			SETVOLTAGE (<i>FuGProbusVRegisterGroups attribute</i>), 119
set_jog_speed() (<i>ILS2T method</i>),	174		SEVENBITS (<i>SerialCommunicationBytesize attribute</i>), 23
set_lower_limits() (<i>PSI9000 method</i>),	108		SHORT_CIRCUIT (<i>NewportSMC100PP.MotorErrors attribute</i>), 146
set_max_acceleration() (<i>ILS2T method</i>),	174		shunt (<i>LEM4000S attribute</i>), 186
set_max_deceleration() (<i>ILS2T method</i>),	174		
set_max_rpm() (<i>ILS2T method</i>),	175		
set_measuring_options() (<i>MBW973 method</i>),	142		
set_message_board() (<i>SupercubeBase method</i>),	40		
set_motor_configuration() (<i>NewportSMC100PP method</i>),			
150			
set_negative_software_limit() (<i>NewportSMC100PP method</i>),	150		
set_number_of_recordings() (<i>HeinzingerDI method</i>),	125		

SHUTDOWN_CURRENT_LIMIT (*PSI9000 attribute*), 107
 SHUTDOWN_VOLTAGE_LIMIT (*PSI9000 attribute*), 107
 ShutterStatus (*CryLasLaser attribute*), 101
 ShutterStatus (*CryLasLaserConfig attribute*), 103
 signal_type (*TiePieGeneratorConfig property*), 89
 SINE (*TiePieGeneratorSignalType attribute*), 89
 SingleCommDevice (*class in hvl_ccb.dev.base*), 97
 SIXBITS (*SerialCommunicationBytesize attribute*), 23
 SIXTEEN (*HeinzingerConfig.RecordingsEnum attribute*), 123
 SIXTEEN_BIT (*TiePieOscilloscopeResolution attribute*), 93
 SL (*NewportConfigCommands attribute*), 145
 SN (*FuGReadbackChannels attribute*), 120
 SOFTWARE_INTERNAL_SIXTY (*CryLasLaser.RepetitionRates attribute*), 101
 SOFTWARE_INTERNAL_TEN (*CryLasLaser.RepetitionRates attribute*), 101
 SOFTWARE_INTERNAL_TWENTY (*CryLasLaser.RepetitionRates attribute*), 101
 software_revision (*LuminoxMeasurementType attribute*), 180
 SPACE (*SerialCommunicationParity attribute*), 25
 SPECIALRAMPUPWARDS (*FuGRampModes attribute*), 120
 spoll() (*VisaCommunication method*), 30
 spoll_handler() (*VisaDevice method*), 184
 SQUARE (*TiePieGeneratorSignalType attribute*), 89
 SR (*NewportConfigCommands attribute*), 145
 srq_mask (*FuGProbusVConfigRegisters property*), 117
 srq_status (*FuGProbusVConfigRegisters property*), 118
 stage_configuration (*NewportSMC100PPConfig attribute*), 153
 START (*LaudaProRp245eCommand attribute*), 137
 start() (*CryLasAttenuator method*), 98
 start() (*CryLasLaser method*), 102
 start() (*Device method*), 95
 start() (*DeviceSequenceMixin method*), 96
 start() (*ExperimentManager method*), 193
 start() (*FuG method*), 112
 start() (*FuGProbusIV method*), 116
 start() (*HeinzingerDI method*), 125
 start() (*HeinzingerPNC method*), 126
 start() (*ILS2T method*), 175
 start() (*LabJack method*), 134
 start() (*LaudaProRp245e method*), 136
 start() (*Luminox method*), 179
 start() (*MBW973 method*), 142
 start() (*NewportSMC100PP method*), 151
 start() (*PfeifferTPG method*), 160
 start() (*PSI9000 method*), 109
 start() (*RTO1024 method*), 170
 start() (*SingleCommDevice method*), 97
 start() (*Supercube2015Base method*), 64
 start() (*SupercubeBase method*), 41
 start() (*Technix method*), 81
 start() (*TiePieGeneratorMixin method*), 89
 start() (*TiePieI2CHostMixin method*), 90
 start() (*TiePieOscilloscope method*), 92
 start() (*VisaDevice method*), 184
 start_control() (*MBW973 method*), 142
 start_measurement() (*TiePieOscilloscope method*), 92
 start_polling() (*Poller method*), 183
 start_ramp() (*LaudaProRp245e method*), 136
 STARTING (*ExperimentStatus attribute*), 193
 States (*NewportSMC100PP attribute*), 146
 status (*ExperimentManager property*), 193
 status (*FuGProbusVConfigRegisters property*), 118
 status (*FuGProbusVDORegisters property*), 118
 status (*Safety attribute*), 56
 status (*Technix property*), 81
 status() (*EarthingStick class method*), 51
 status_1 (*Door attribute*), 49
 status_1 (*EarthingRod attribute*), 49
 status_1 (*EarthingStick attribute*), 51
 status_1_closed (*EarthingStick attribute*), 68
 status_1_connected (*EarthingStick attribute*), 68
 status_1_open (*EarthingStick attribute*), 68
 status_2 (*Door attribute*), 49
 status_2 (*EarthingRod attribute*), 49
 status_2 (*EarthingStick attribute*), 51
 status_2_closed (*EarthingStick attribute*), 69
 status_2_connected (*EarthingStick attribute*), 69
 status_2_open (*EarthingStick attribute*), 69
 status_3 (*Door attribute*), 49
 status_3 (*EarthingRod attribute*), 50
 status_3 (*EarthingStick attribute*), 51
 status_3_closed (*EarthingStick attribute*), 69
 status_3_connected (*EarthingStick attribute*), 69
 status_3_open (*EarthingStick attribute*), 69
 status_4 (*EarthingStick attribute*), 51
 status_4_closed (*EarthingStick attribute*), 69
 status_4_connected (*EarthingStick attribute*), 69
 status_4_open (*EarthingStick attribute*), 69
 status_5 (*EarthingStick attribute*), 51
 status_5_closed (*EarthingStick attribute*), 69
 status_5_connected (*EarthingStick attribute*), 69
 status_5_open (*EarthingStick attribute*), 69
 status_6 (*EarthingStick attribute*), 51
 status_6_closed (*EarthingStick attribute*), 69
 status_6_connected (*EarthingStick attribute*), 69
 status_6_open (*EarthingStick attribute*), 69
 status_closed() (*EarthingStick class method*), 69
 status_connected() (*EarthingStick class method*), 69
 status_error (*Safety attribute*), 73
 status_green (*Safety attribute*), 73
 status_open() (*EarthingStick class method*), 69

status_ready_for_red (*Safety attribute*), 73
status_red (*Safety attribute*), 73
STATUSBYTE (*FuGReadbackChannels attribute*), 120
statuses() (*EarthingStick class method*), 51
stop (*Errors attribute*), 52, 70
STOP (*LaudaProRp245eCommand attribute*), 137
stop() (*CryLasLaser method*), 102
stop() (*Device method*), 95
stop() (*DeviceSequenceMixin method*), 96
stop() (*ExperimentManager method*), 193
stop() (*FuGProbusIV method*), 116
stop() (*HeinzingerDI method*), 125
stop() (*ILS2T method*), 175
stop() (*LabJack method*), 134
stop() (*LaudaProRp245e method*), 136
stop() (*Luminox method*), 179
stop() (*MBW973 method*), 142
stop() (*NewportSMC100PP method*), 151
stop() (*PfeifferTPG method*), 161
stop() (*PSI9000 method*), 109
stop() (*RTO1024 method*), 170
stop() (*SingleCommDevice method*), 97
stop() (*Supercube2015Base method*), 64
stop() (*SupercubeBase method*), 41
stop() (*Technix method*), 81
stop() (*TiePieGeneratorMixin method*), 89
stop() (*TiePieI2CHostMixin method*), 90
stop() (*TiePieOscilloscope method*), 92
stop() (*VisaDevice method*), 184
stop_acquisition() (*RTO1024 method*), 170
stop_motion() (*NewportSMC100PP method*), 151
stop_number (*Errors attribute*), 70
stop_polling() (*Poller method*), 183
stop_ramp() (*LaudaProRp245e method*), 136
stopbits (*CryLasAttenuatorSerialCommunicationConfig attribute*), 100
stopbits (*CryLasLaserSerialCommunicationConfig attribute*), 106
stopbits (*FuGSerialCommunicationConfig attribute*), 122
stopbits (*HeinzingerSerialCommunicationConfig attribute*), 127
stopbits (*LuminoxSerialCommunicationConfig attribute*), 182
stopbits (*MBW973SerialCommunicationConfig attribute*), 144
stopbits (*NewportSMC100PPSerialCommunicationConfig attribute*), 157
stopbits (*PfeifferTPGSerialCommunicationConfig attribute*), 163
Stopbits (*SerialCommunicationConfig attribute*), 24
stopbits (*SerialCommunicationConfig attribute*), 25
streaming (*LuminoxOutputMode attribute*), 181
StrEnumBase (*class in hvl_ccb.utils.enum*), 188
sub_handler (*OpcUaCommunicationConfig attribute*), 22
sub_handler (*SupercubeOpcUaCommunicationConfig attribute*), 43, 66
suitable_range() (*TiePieOscilloscopeRange static method*), 86
Supercube2015Base (*class in hvl_ccb.dev.supercube2015.base*), 62
Supercube2015WithFU (*class in hvl_ccb.dev.supercube2015.typ_a*), 74
SupercubeA0pcUaCommunication (*class in hvl_ccb.dev.supercube.typ_a*), 57
SupercubeA0pcUaCommunication (*class in hvl_ccb.dev.supercube2015.typ_a*), 75
SupercubeA0pcUaConfiguration (*class in hvl_ccb.dev.supercube.typ_a*), 57
SupercubeA0pcUaConfiguration (*class in hvl_ccb.dev.supercube2015.typ_a*), 75
SupercubeB (*class in hvl_ccb.dev.supercube.typ_b*), 60
SupercubeBase (*class in hvl_ccb.dev.supercube.base*), 37
SupercubeB0pcUaCommunication (*class in hvl_ccb.dev.supercube.typ_b*), 60
SupercubeB0pcUaConfiguration (*class in hvl_ccb.dev.supercube.typ_b*), 60
SupercubeConfiguration (*class in hvl_ccb.dev.supercube.base*), 41
SupercubeConfiguration (*class in hvl_ccb.dev.supercube2015.base*), 65
SupercubeEarthlingStickOperationError, 42
SupercubeOpcEndpoint (*class in hvl_ccb.dev.supercube.constants*), 56
SupercubeOpcEndpoint (*class in hvl_ccb.dev.supercube2015.constants*), 74
SupercubeOpcUaCommunication (*class in hvl_ccb.dev.supercube.base*), 42
SupercubeOpcUaCommunication (*class in hvl_ccb.dev.supercube2015.base*), 65
SupercubeOpcUaCommunicationConfig (*class in hvl_ccb.dev.supercube.base*), 42
SupercubeOpcUaCommunicationConfig (*class in hvl_ccb.dev.supercube2015.base*), 65
SupercubeSubscriptionHandler (*class in hvl_ccb.dev.supercube.base*), 43
SupercubeSubscriptionHandler (*class in hvl_ccb.dev.supercube2015.base*), 66
SupercubeWithFU (*class in hvl_ccb.dev.supercube.typ_a*), 58
switch_to_operate (*Safety attribute*), 56
switch_to_ready (*Safety attribute*), 56
switchto_green (*Safety attribute*), 73
switchto_operate (*Safety attribute*), 73
switchto_ready (*Safety attribute*), 73
SyncCommunicationProtocol (*class in*

hvl_ccb.comm.base), 14
SyncCommunicationProtocolConfig (class in
hvl_ccb.comm.base), 14

T

T (*LabJack.ThermocoupleType attribute), 131*
t13_1 (*GeneralSockets attribute), 52, 70*
t13_2 (*GeneralSockets attribute), 52, 70*
t13_3 (*GeneralSockets attribute), 52, 70*
T13_SOCKET_PORTS (in module
hvl_ccb.dev.supercube.constants), 57
T13_SOCKET_PORTS (in module
hvl_ccb.dev.supercube2015.constants), 74
T1MS (*FuGMonitorModes attribute), 115*
T200MS (*FuGMonitorModes attribute), 115*
T20MS (*FuGMonitorModes attribute), 115*
T256US (*FuGMonitorModes attribute), 115*
T4 (*LabJack.DeviceType attribute), 130*
T4 (*LJMCommunicationConfig.DeviceType attribute), 16
T40MS (*FuGMonitorModes attribute), 115*
T4MS (*FuGMonitorModes attribute), 115*
T560 (class in *hvl_ccb.dev.higland_t560.device*), 35
T560Communication (class in
hvl_ccb.dev.higland_t560.base), 33
T560CommunicationConfig (class in
hvl_ccb.dev.higland_t560.base), 33
T560Config (class in *hvl_ccb.dev.higland_t560.device*),
36
T560Error, 34
T7 (*LabJack.DeviceType attribute), 130*
T7 (*LJMCommunicationConfig.DeviceType attribute), 16
T7_PRO (*LabJack.DeviceType attribute), 130*
T7_PRO (*LJMCommunicationConfig.DeviceType attribute), 16
T800MS (*FuGMonitorModes attribute), 115*
T80MS (*FuGMonitorModes attribute), 116*
target_pulse_energy (*CryLasLaser property), 103*
Tcp (class in *hvl_ccb.comm.tcp*), 26
TCP (*LJMCommunicationConfig.ConnectionType attribute), 16
TcpCommunicationConfig (class in *hvl_ccb.comm.tcp*),
26
TCPIP_INSTR (*VisaCommunicationConfig.InterfaceType attribute), 30*
TCPIP_SOCKET (*VisaCommunicationConfig.InterfaceType attribute), 30*
TEC1 (*CryLasLaser.AnswersStatus attribute), 101*
TEC2 (*CryLasLaser.AnswersStatus attribute), 101*
Technix (class in *hvl_ccb.dev.technix.device*), 80
TechnixConfig (class in *hvl_ccb.dev.technix.device*), 81
TechnixError, 77
TechnixFaultError, 77
TechnixSerialCommunication (class in
hvl_ccb.dev.technix.base), 77
TechnixSerialCommunicationConfig (class in
hvl_ccb.dev.technix.base), 77
TechnixTelnetCommunication (class in
hvl_ccb.dev.technix.base), 78
TechnixTelnetCommunicationConfig (class in
hvl_ccb.dev.technix.base), 78
TelnetCommunication (class in *hvl_ccb.comm.telnet*),
27
TelnetCommunicationConfig (class in
hvl_ccb.comm.telnet), 28
TelnetError, 29
TEMP (*ILS2TRegAddr attribute), 177*
TEMP_SET_POINT (*LaudaProRp245eCommand attribute), 137
temp_set_point_init (*LaudaProRp245eConfig attribute), 139
Temperature (class in *hvl_ccb.utils.conversion_unit*),
187
temperature_sensor (*LuminoxMeasurementType attribute), 181
temperature_unit (*LMT70A attribute), 186
TEN (*LabJack.AInRange attribute), 129*
TEN (*LabJack.CalMicroAmpere attribute), 129*
TEN_MHZ (*LabJack.ClockFrequency attribute), 129*
terminator (*AsyncCommunicationProtocolConfig attribute), 14
terminator (*CryLasAttenuatorSerialCommunicationConfig attribute), 100
terminator (*CryLasLaserSerialCommunicationConfig attribute), 106
TERMINATOR (*FuGProbusIVCommands attribute), 117
terminator (*FuGProbusVConfigRegisters property),
118
terminator (*FuGSerialCommunicationConfig attribute), 122
terminator (*HeinzingerSerialCommunicationConfig attribute), 127
terminator (*LaudaProRp245eTcpCommunicationConfig attribute), 140
terminator (*LuminoxSerialCommunicationConfig attribute), 182
terminator (*MBW973SerialCommunicationConfig attribute), 144
terminator (*NewportSMC100PPSSerialCommunicationConfig attribute), 157
terminator (*PfeifferTPGSerialCommunicationConfig attribute), 163
terminator (*T560CommunicationConfig attribute), 34
terminator_str() (*SerialCommunicationConfig method), 25
THIRTY_TWO_BIT (*LabJack.BitLimit attribute), 129
TiePieDeviceConfig (class in *hvl_ccb.dev.tiepie.base*),
83
TiePieDeviceType (class in *hvl_ccb.dev.tiepie.base*),***********************

84		
TiePieError, 84		
TiePieGeneratorConfig (class <i>hvl_ccb.dev.tiepie.generator</i>), 88	in	attribute), 157
TiePieGeneratorConfigLimits (class <i>hvl_ccb.dev.tiepie.generator</i>), 89	in	timeout (<i>PfeifferTPGSerialCommunicationConfig</i> <i>attribute</i>), 163
TiePieGeneratorMixin (class <i>hvl_ccb.dev.tiepie.generator</i>), 89	in	timeout (<i>SerialCommunicationConfig</i> <i>attribute</i>), 25
TiePieGeneratorSignalType (class <i>hvl_ccb.dev.tiepie.generator</i>), 89	in	timeout (<i>TelnetCommunicationConfig</i> <i>attribute</i>), 29
TiePieHS5 (class in <i>hvl_ccb.dev.tiepie.device</i>), 87		timeout (<i>VisaCommunicationConfig</i> <i>attribute</i>), 32
TiePieHS6 (class in <i>hvl_ccb.dev.tiepie.device</i>), 88		Torr (<i>PfeifferTPG.PressureUnits</i> <i>attribute</i>), 158
TiePieI2CHostConfig (class in <i>hvl_ccb.dev.tiepie.i2c</i>), 90		TORR (<i>Pressure</i> <i>attribute</i>), 187
TiePieI2CHostConfigLimits (class <i>hvl_ccb.dev.tiepie.i2c</i>), 90	in	TPG25xA (<i>PfeifferTPGConfig.Model</i> <i>attribute</i>), 161
TiePieI2CHostMixin (class in <i>hvl_ccb.dev.tiepie.i2c</i>), 90		TPGx6x (<i>PfeifferTPGConfig.Model</i> <i>attribute</i>), 161
TiePieOscilloscope (class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 91	in	TPR (<i>PfeifferTPG.SensorTypes</i> <i>attribute</i>), 159
TiePieOscilloscopeAutoResolutionModes (class in <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 92		transmission (<i>CryLasAttenuator</i> <i>property</i>), 98
TiePieOscilloscopeChannelConfig (class <i>hvl_ccb.dev.tiepie.channel</i>), 85	in	TRIANGLE (<i>TiePieGeneratorSignalType</i> <i>attribute</i>), 89
TiePieOscilloscopeChannelConfigLimits (class in <i>hvl_ccb.dev.tiepie.channel</i>), 86		trigger_enabled (<i>TiePieOscilloscopeChannelConfig</i> <i>property</i>), 86
TiePieOscilloscopeChannelCoupling (class <i>hvl_ccb.dev.tiepie.channel</i>), 86	in	trigger_hysteresis (<i>TiePieOscilloscopeChannel-</i> <i>Config</i> <i>property</i>), 86
TiePieOscilloscopeConfig (class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 92	in	trigger_kind (<i>TiePieOscilloscopeChannelConfig</i> <i>prop-</i> <i>erty</i>), 86
TiePieOscilloscopeConfigLimits (class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 93	in	trigger_level (<i>T560</i> <i>property</i>), 36
TiePieOscilloscopeRange (class <i>hvl_ccb.dev.tiepie.channel</i>), 86	in	trigger_level (<i>TiePieOscilloscopeChannelConfig</i> <i>property</i>), 86
TiePieOscilloscopeResolution (class <i>hvl_ccb.dev.tiepie.oscilloscope</i>), 93	in	trigger_level_mode (<i>TiePieOscilloscopeChannel-</i> <i>Config</i> <i>property</i>), 86
TiePieOscilloscopeTriggerKind (class <i>hvl_ccb.dev.tiepie.channel</i>), 86	in	trigger_mode (<i>T560</i> <i>property</i>), 36
TiePieOscilloscopeTriggerLevelMode (class <i>hvl_ccb.dev.tiepie.channel</i>), 87	in	trigger_timeout (<i>TiePieOscilloscopeConfig</i> <i>prop-</i> <i>erty</i>), 93
TiePieWS5 (class in <i>hvl_ccb.dev.tiepie.device</i>), 88		triggered (<i>BreakdownDetection</i> <i>attribute</i>), 49, 68
timeout (<i>CryLasAttenuatorSerialCommunicationConfig</i> <i>attribute</i>), 100		TriggerMode (class in <i>hvl_ccb.dev.highland_t560.base</i>), 34
timeout (<i>CryLasLaserSerialCommunicationConfig</i> <i>at-</i> <i>tribute</i>), 106		TWELVE_BIT (<i>TiePieOscilloscopeResolution</i> <i>attribute</i>), 93
timeout (<i>FuGSerialCommunicationConfig</i> <i>attribute</i>), 122		TWELVE_HUNDRED_FIFTY_KHZ (<i>Lab-</i> <i>Jack.ClockFrequency</i> <i>attribute</i>), 129
timeout (<i>HeinzingerSerialCommunicationConfig</i> <i>attribute</i>), 127		TWENTY_FIVE_HUNDRED_KHZ (<i>LabJack.ClockFrequency</i> <i>attribute</i>), 129
timeout (<i>LuminoxSerialCommunicationConfig</i> <i>at-</i> <i>tribute</i>), 182		TWENTY_MHZ (<i>LabJack.ClockFrequency</i> <i>attribute</i>), 129
timeout (<i>MBW973SerialCommunicationConfig</i> <i>at-</i> <i>tribute</i>), 144		TWENTY_VOLT (<i>TiePieOscilloscopeRange</i> <i>attribute</i>), 86
timeout (<i>NewportSMC100PPSerialCommunicationConfig</i>		U
		Underrange (<i>PfeifferTPG.SensorStatus</i> <i>attribute</i>), 159
		Unit (class in <i>hvl_ccb.utils.conversion_unit</i>), 187
		unit (<i>ILS2TModbusTcpCommunicationConfig</i> <i>attribute</i>), 177
		unit (<i>ModbusTcpCommunicationConfig</i> <i>attribute</i>), 19
		unit (<i>PfeifferTPG</i> <i>property</i>), 161
		unit_current (<i>HeinzingerPNC</i> <i>property</i>), 126
		unit_voltage (<i>HeinzingerPNC</i> <i>property</i>), 126

UNKNOWN (*HeinzingerDI.OutputStatus attribute*), 124
 UNKNOWN (*HeinzingerPNC.UnitCurrent attribute*), 125
 UNKNOWN (*HeinzingerPNC.UnitVoltage attribute*), 125
 UNKNOWN (*TiePieGeneratorSignalType attribute*), 89
 UNKNOWN (*TiePieOscilloscopeAutoResolutionModes attribute*), 92
 UNKNOWN (*TiePieOscilloscopeTriggerLevelMode attribute*), 87
 UNREADY_INACTIVE (*CryLasLaser.LaserStatus attribute*), 101
 update_laser_status() (*CryLasLaser method*), 103
 update_period (*OpcUaCommunicationConfig attribute*), 22
 update_repetition_rate() (*CryLasLaser method*), 103
 update_shutter_status() (*CryLasLaser method*), 103
 update_target_pulse_energy() (*CryLasLaser method*), 103
 UpdateEspStageInfo (*NewportSMC100PPConfig.EspStageConfig attribute*), 152
 UPPER_TEMP (*LaudaProRp245eCommand attribute*), 137
 upper_temp (*LaudaProRp245eConfig attribute*), 139
 USB (*LaudaProRp245eConfig.ExtControlModeEnum attribute*), 138
 USB (*LJMCommunicationConfig.ConnectionType attribute*), 16
 use_external_clock() (*T560 method*), 36
 user_position_offset (*NewportSMC100PPConfig attribute*), 153
 user_steps() (*ILS2T method*), 175

V

V (*HeinzingerPNC.UnitVoltage attribute*), 125
 VA (*NewportConfigCommands attribute*), 145
 validate_and_resolve_host() (*in module hvl_ccb.utils.validation*), 189
 validate_bool() (*in module hvl_ccb.utils.validation*), 189
 validate_number() (*in module hvl_ccb.utils.validation*), 189
 validate_pump_level() (*LaudaProRp245e method*), 136
 validate_tcp_port() (*in module hvl_ccb.utils.validation*), 190
 value (*FuGProbusVMonitorRegisters property*), 119
 value (*LabJack.AInRange property*), 129
 value_raw (*FuGProbusVMonitorRegisters property*), 119
 ValueEnum (*class in hvl_ccb.utils.enum*), 188
 VB (*NewportConfigCommands attribute*), 145
 velocity (*NewportSMC100PPConfig attribute*), 153
 visa_backend (*VisaCommunicationConfig attribute*), 32

VisaCommunication (*class in hvl_ccb.comm.visa*), 29
 VisaCommunicationConfig (*class in hvl_ccb.comm.visa*), 30
 VisaCommunicationConfig.InterfaceType (*class in hvl_ccb.comm.visa*), 30
 VisaCommunicationError, 32
 VisaDevice (*class in hvl_ccb.dev.visa*), 183
 VisaDeviceConfig (*class in hvl_ccb.dev.visa*), 184
 VOLT (*ILS2TRegAddr attribute*), 177
 Volt (*PfeifferTPG.PressureUnits attribute*), 158
 voltage (*FuG property*), 113
 VOLTAGE (*FuGProbusIVCommands attribute*), 117
 VOLTAGE (*FuGReadbackChannels attribute*), 120
 voltage (*Technix property*), 81
 voltage_lower_limit (*PSI9000Config attribute*), 110
 voltage_max (*Power attribute*), 55, 72
 voltage_monitor (*FuG property*), 113
 voltage_primary (*Power attribute*), 55, 72
 voltage_regulation (*Technix property*), 81
 voltage_slope (*Power attribute*), 55, 72
 voltage_target (*Power attribute*), 55, 72
 voltage_upper_limit (*PSI9000Config attribute*), 110

W

WAIT_AFTER_WRITE (*VisaCommunication attribute*), 30
 wait_for_polling_result() (*Poller method*), 183
 wait_operation_complete() (*VisaDevice method*), 184
 wait_sec_initialisation (*PSI9000Config attribute*), 110
 wait_sec_max_disable (*ILS2TConfig attribute*), 176
 wait_sec_post_absolute_position (*ILS2TConfig attribute*), 176
 wait_sec_post_activate (*LuminoxConfig attribute*), 180
 wait_sec_post_cannot_disable (*ILS2TConfig attribute*), 176
 wait_sec_post_enable (*ILS2TConfig attribute*), 176
 wait_sec_post_relative_step (*ILS2TConfig attribute*), 176
 wait_sec_pre_read_or_write (*LaudaProRp245eTcpCommunicationConfig attribute*), 140
 wait_sec_read_text_nonempty (*AsyncCommunicationProtocolConfig attribute*), 14
 wait_sec_read_text_nonempty (*FuGSerialCommunicationConfig attribute*), 122
 wait_sec_read_text_nonempty (*HeinzingerSerialCommunicationConfig attribute*), 128
 wait_sec_retry_get_device (*TiePieDeviceConfig attribute*), 84
 wait_sec_settings_effect (*PSI9000Config attribute*), 110
 wait_sec_stop_commands (*FuGConfig attribute*), 113

wait_sec_stop_commands (*HeinzingerConfig attribute*), 123
wait_sec_system_lock (*PSI9000Config attribute*), 110
wait_sec_trials_activate (*LuminoxConfig attribute*), 180
wait_timeout_retry_sec (*OpcUaCommunicationConfig attribute*), 22
wait_until_motor_initialized() (*NewportSMC100PP method*), 151
wait_until_ready() (*CryLasLaser method*), 103
warning (*Errors attribute*), 52
waveform (*TiePieGeneratorConfig property*), 89
WIFI (*LJMCommunicationConfig.ConnectionType attribute*), 16
wrap_libtiepie_exception() (*in module hvl_ccb.dev.tiepie.base*), 84
write() (*AsyncCommunicationProtocol method*), 12
write() (*MBW973 method*), 142
write() (*OpcUaCommunication method*), 21
write() (*Supercube2015Base method*), 64
write() (*SupercubeBase method*), 41
write() (*Tcp method*), 26
write() (*VisaCommunication method*), 30
write_absolute_position() (*ILS2T method*), 175
write_bytes() (*AsyncCommunicationProtocol method*), 12
write_bytes() (*SerialCommunication method*), 23
write_bytes() (*TelnetCommunication method*), 28
write_command() (*LauraProRp245eTcpCommunication method*), 140
write_name() (*LJMCommunication method*), 16
write_names() (*LJMCommunication method*), 16
write_registers() (*ModbusTcpCommunication method*), 18
write_relative_step() (*ILS2T method*), 175
write_termination (*VisaCommunicationConfig attribute*), 32
write_text() (*AsyncCommunicationProtocol method*), 13
WRONG_ESP_STAGE (*NewportSMC100PP.MotorErrors attribute*), 146

X

x_stat (*FuGProbusVDIRegisters property*), 118
XOUTPUTS (*FuGProbusIVCommands attribute*), 117

Y

YES (*FuGDigitalVal attribute*), 114

Z

ZX (*NewportConfigCommands attribute*), 145